



Miltenyi Biotec

CliniMACS[®] Cell Separation System

Product catalog 2008

Product overview

Page Product Regulatory status¹⁾ Availability Order no.

CliniMACS® Plus Instrument

7	CliniMACS Plus Instrument	CE	Europe, CAN, AUS, US	151-01
7	CliniMACS Plus Upgrade Kit "Software"		Europe, CAN, AUS, US ²⁾	755-16

CliniMACS® Service Contracts

8	CliniMACS Full Service Agreement		Europe	160-000-973
8	CliniMACS Maintenance		Europe	160-000-974
8	CliniMACS Premium Service Agreement		CAN, US	160-000-701
8	CliniMACS Basic Service Agreement		CAN, US	160-000-702

CliniMACS® Tubing Sets

9	CliniMACS Tubing Set	CE	Europe, CAN, AUS, US ²⁾	161-01
9	CliniMACS Tubing Set LS	CE	Europe, CAN, AUS, US ²⁾	162-01
10	CliniMACS Tubing Set for Research Use	For research use only, not for use in humans	US	165-01
10	CliniMACS Tubing Set LS for Research Use	For research use only, not for use in humans	US	168-01
11	CliniMACS Depletion Tubing Set	CE	Europe, US ²⁾	261-01
11	CliniMACS Depletion Tubing Set for Research Use	For research use only, not for use in humans	US, CAN, AUS	266-01
12	CliniMACS Tubing Set 600 for Research Use	For research use only, not for use in humans	Worldwide	166-01
12	CliniMACS Tubing Set 150 for Research Use	For research use only, not for use in humans	Worldwide	167-01

For availability in your country please contact your local representative

1) The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE). CliniMACS MicroBeads are for research use only, and not for use in humans.

2) In the USA, the CliniMACS® System components including the CliniMACS Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

Product overview

Page Product Regulatory status¹⁾ Availability Order no.

Stem and progenitor cells

13	CliniMACS CD34 Reagent	CE	Europe, CAN, AUS, US ²⁾	171-01
13	CliniMACS CD34 Complete Kit	CE-marked components	Europe	178-01
13	CliniMACS CD34 Complete Kit with Accessories	CE-marked components	Europe	177-01
13	CliniMACS CD34 Complete LS Kit	CE-marked components	Europe	170-01
13	CliniMACS CD34 Complete LS Kit with Accessories	CE-marked components	Europe	169-01
13	CliniMACS CD34 MicroBeads	For research use only, not for use in humans	US	175-01
16	CliniMACS CD133 Reagent	CE	Europe, US ²⁾	172-01
16	CliniMACS CD133 Complete Kit	CE-marked components	Europe	196-01
16	CliniMACS CD133 Complete Kit with Accessories	CE-marked components	Europe	198-01
16	CliniMACS CD133 Complete LS Kit	CE-marked components	Europe	197-01
16	CliniMACS CD133 Complete LS Kit with Accessories	CE-marked components	Europe	199-01
16	CliniMACS CD133 MicroBeads	For research use only, not for use in humans	US, CAN, AUS	195-01
18	CliniMACS CD3/CD19 Reagent Combination	CE-marked components	Europe	704-00
18	CliniMACS CD3/CD19 Complete Kit	CE-marked components	Europe	740-01
18	CliniMACS CD3/CD19 Large-Scale Complete Kit	CE-marked components	Europe	740-02

T cells

20	CliniMACS CD3 Reagent	CE	Europe, US ²⁾	273-01
20	CliniMACS CD3 MicroBeads	For research use only, not for use in humans	US, CAN, AUS	176-01
21	CliniMACS CD4 Reagent	CE	Europe, US ²⁾	276-01
21	CliniMACS CD4 MicroBeads	For research use only, not for use in humans	US, CAN, AUS	304-01
22	CliniMACS CD8 Reagent	CE	Europe, US ²⁾	275-01
22	CliniMACS CD8 MicroBeads	For research use only, not for use in humans	US, CAN, AUS	308-01
23	CliniMACS CD25 Reagent	CE	Europe, US ²⁾	NEW 274-01
23	CliniMACS CD25 MicroBeads	For research use only, not for use in humans	US, CAN, AUS	325-01

CD3/CD19 combination products, please see chapter "Stem and progenitor cells"

Antigen-specific T cells

25	CliniMACS Cytokine Capture System (IFN-gamma)	CE	Europe, US ²⁾	279-01
25	CliniMACS Cytokine Capture System (IFN-gamma) MicroBeads	For research use only, not for use in humans	US, CAN, AUS	287-01

For availability in your country please contact your local representative

1) The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE). CliniMACS MicroBeads are for research use only, and not for use in humans.

2) In the USA, the CliniMACS® System components including the CliniMACS Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

Product overview

Page	Product	Regulatory status ¹⁾	Availability	Order no.
NK cells				
28	CliniMACS CD56 Reagent	CE	Europe, US ²⁾	271-01
28	CliniMACS CD3/CD56 Reagent Combination	CE-marked components	Europe	704-01
28	CliniMACS CD3/CD56 Complete Kit	CE-marked components	Europe	740-04
28	CliniMACS CD56 MicroBeads	For research use only, not for use in humans	US, CAN, AUS	194-01
B cells				
30	CliniMACS CD19 Reagent	CE	Europe, US ²⁾	179-01
30	CliniMACS CD19 MicroBeads	For research use only, not for use in humans	US, CAN, AUS	193-01
Dendritic cells				
32	CliniMACS CD14 Reagent	CE	Europe, US ²⁾	272-01
32	CliniMACS CD14 Complete Kit	CE-marked components	Europe	281-01
32	CliniMACS CD14 MicroBeads	For research use only, not for use in humans	US, CAN, AUS	191-01
33	CliniMACS CD1c (BDCA-1)-Biotin	CE	Europe, US ²⁾	NEW 277-01
33	CliniMACS CD1c (BDCA-1)/CD19 Complete Kit	CE-marked component	Europe	NEW 740-03
33	CliniMACS CD1c (BDCA-1)-Biotin for Research Use	For research use only, not for use in humans	US, CAN, AUS	255-01
34	CliniMACS CD304 (BDCA-4) MicroBeads	For research use only, not for use in humans	Worldwide	291-01
Flexible labeling system				
35	CliniMACS Anti-Biotin Reagent	CE	Europe, US ²⁾	173-01
35	CliniMACS Anti-Biotin MicroBeads	For research use only, not for use in humans	US, CAN, AUS	192-01
35	CliniMACS® Biotinylation Kit 500	CE	Europe, US ²⁾	NEW 701-06
35	CliniMACS® Biotinylation Kit 500 for Research Use	For research use only, not for use in humans	US, CAN, AUS	257-01
CliniMACS® Buffers				
36	CliniMACS PBS/EDTA Buffer	CE	Europe, CAN, AUS, US ²⁾	700-25
36	CliniMACS PBS/EDTA Buffer for Research Use	For research use only, not for use in humans	US	705-25

For availability in your country please contact your local representative

1) The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE). CliniMACS MicroBeads are for research use only, and not for use in humans.

2) In the USA, the CliniMACS® System components including the CliniMACS Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

Product overview

Page Product Regulatory status¹⁾ Availability Order no.

Accessories

38	Cell Expansion Bag (tube)	CE ³⁾	Europe	200-074-301
38	Cell Expansion Bag (6 ports)	CE ³⁾	Europe	200-074-302
38	Cell Expansion Bag (tube)	For research use only	US, CAN, AUS	NEW 130-074-351
38	Cell Expansion Bag (6 ports)	For research use only	US, CAN, AUS	NEW 130-074-352
39	Cell Differentiation Bag (tube) – 100	CE ³⁾	Europe	200-074-101
39	Cell Differentiation Bag (tube) – 250	CE ³⁾	Europe	200-074-102
39	Cell Differentiation Bag (tube) – 500	CE ³⁾	Europe	200-074-103
39	Cell Differentiation Bag (6 ports) – 100	CE ³⁾	Europe	200-074-201
39	Cell Differentiation Bag (6 ports) – 250	CE ³⁾	Europe	200-074-202
39	Cell Differentiation Bag (6 ports) – 500	CE ³⁾	Europe	200-074-203
39	Cell Differentiation Bag (tube) – 100	For research use only	US, CAN, AUS	NEW 130-074-151
39	Cell Differentiation Bag (tube) – 250	For research use only	US, CAN, AUS	NEW 130-074-152
39	Cell Differentiation Bag (tube) – 500	For research use only	US, CAN, AUS	NEW 130-074-153
39	Cell Differentiation Bag (6 ports) – 100	For research use only	US, CAN, AUS	NEW 130-074-251
39	Cell Differentiation Bag (6 ports) – 250	For research use only	US, CAN, AUS	NEW 130-074-252
39	Cell Differentiation Bag (6 ports) – 500	For research use only	US, CAN, AUS	NEW 130-074-253
40	Pre-System Filter	CE	Europe	181-01
40	Luer/Spike Interconnector	CE	Europe	187-01
40	Sampling Site Coupler	CE	Europe	189-01
41	Transfer Set Coupler/Needle	CE	Europe	185-01
41	Transfer Set Coupler/Coupler	CE	Europe	186-01
41	Transfer Set Needle/Needle	CE	Europe	188-01
42	Transfer Bag 150 mL	CE	Europe	183-01
42	Transfer Bag 600 mL	CE	Europe	190-01
42	Transfer Bag 1000 mL	CE	Europe	180-01
42	Transfer Bag 600 mL with 8 couplers	CE	Europe	184-01

For availability in your country please contact your local representative

1) The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE). CliniMACS MicroBeads are for research use only, and not for use in humans.

2) In the USA, the CliniMACS® System components including the CliniMACS Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

3) The Cell Expansion Bags and the Cell Differentiation Bags are manufactured and controlled under an ISO 13485 certified quality system. These products are available in Europe as CE-marked medical devices.

CliniMACS® Technology

The CliniMACS® Plus Instrument is an automated cell separation system based on MACS® Technology. It enables the operator to perform clinical-scale magnetic enrichment of target cells or depletion of unwanted cells in a closed and sterile system.

The cells are first magnetically labeled in the Cell Preparation Bag using an antigen-specific reagent. After labeling, the cells are washed to remove excess reagent. Then the Cell Preparation Bag containing the labeled cells is connected to the Tubing Set, which in turn is connected to the CliniMACS Buffer Bag and Cell Collection Bag. After starting, for example the enrichment program, the system automatically applies the cell sample onto the selection column. The sample application is followed by a series of washing steps depending on the chosen enrichment program. Finally, the target cells are eluted from the column. The CliniMACS Plus Instrument in combination with the CliniMACS® System separates large numbers of cells providing target cells with high purity and excellent yield. A range of tubing sets optimized for different applications is available.

The CliniMACS System (CliniMACS Plus Instrument, CliniMACS Tubing Sets, CliniMACS Reagents, and CliniMACS PBS/EDTA Buffer) offers a flexible platform enabling the separation of many cell types as well as customized separation protocols according to specific sample requirements. Unique reagents and tubing sets also allow for the transition of promising research applications in innovative cell therapy concepts.



Figure 1: Cell Selection Column placed in the CliniMACS Magnetic Separation Unit.

The CliniMACS Plus Instrument is CE-marked for clinical use in Europe and is used for the separation of a variety of human cell types in combination with a CliniMACS Tubing Set, a CliniMACS Reagent, and the CliniMACS PBS/EDTA Buffer. The portfolio of CliniMACS Reagents includes reagents for the separation

of stem cells, monocytes, DCs, NK cells, B cells, and T cell subsets. In the USA, the CliniMACS® System components including the CliniMACS Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).



Figure 2: The CliniMACS Plus Instrument with installed CliniMACS Tubing Set.

CliniMACS® Plus Instrument


CliniMACS® Plus Instrument

The key components of the CliniMACS® Plus Instrument are the integrated micro-computer, the magnetic separation unit, the peristaltic pump, and various pinch valves. The integrated computer

controls all components of the instrument and directs the system to perform reproducible procedures in a standardized sequence. The magnetic separation unit includes a movable permanent magnet and a holder for the

selection column. The peristaltic pump controls constant flow rates throughout the tubing set and ensures, together with the pinch valves, the controlled flow of buffer and cell suspension.

Key components of the CliniMACS Plus Instrument with installed CliniMACS® Tubing Set.

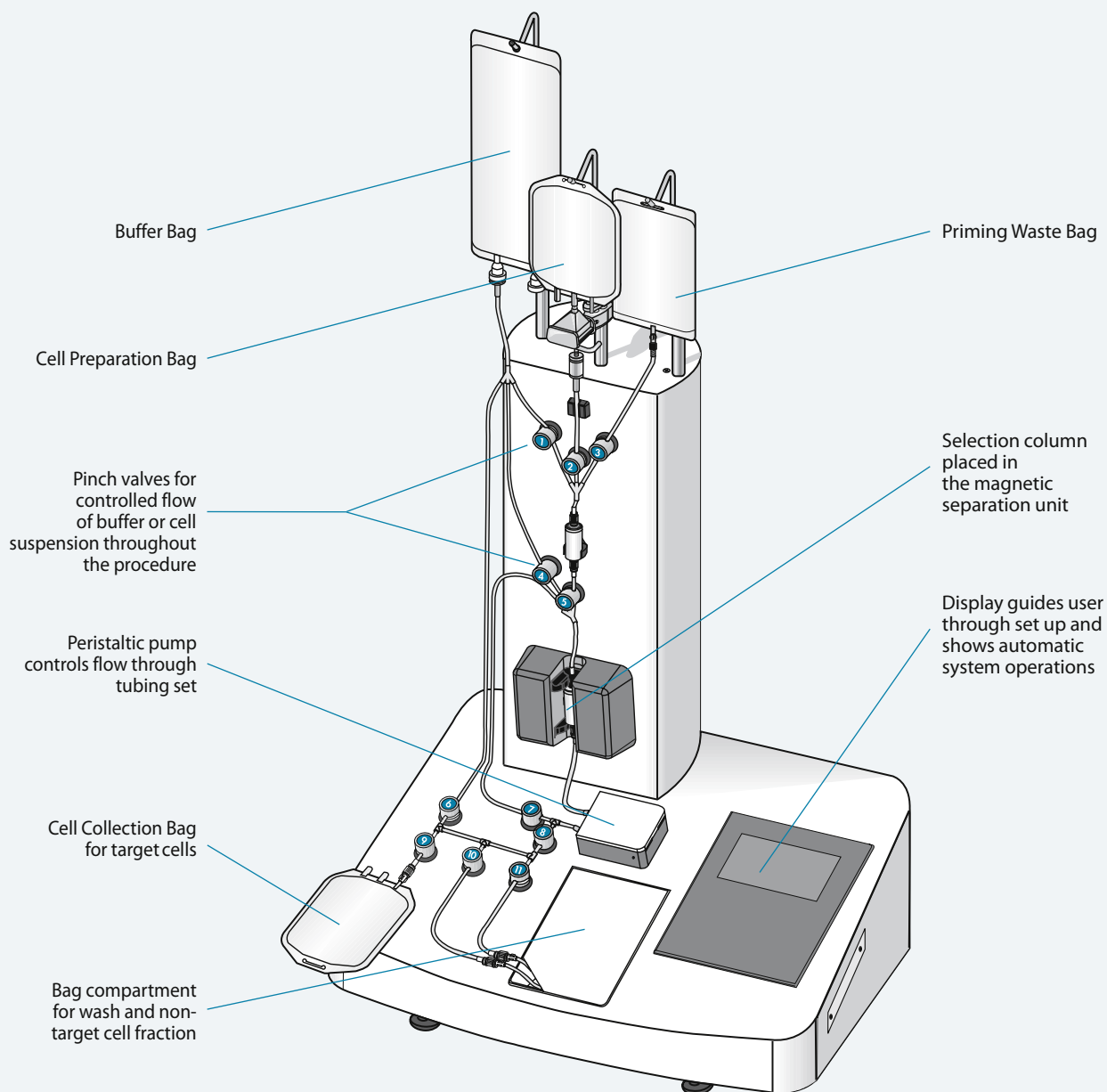

CliniMACS® Plus Instrument

CliniMACS® Plus Instrument			
Product	Regulatory status	Availability	Order no.
CliniMACS Plus Instrument	CE	Europe, CAN, AUS, US ¹⁾	151-01
For availability in your country please contact your local representative.			
1) In the USA, the CliniMACS System components including the CliniMACS Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).			

Description

The CliniMACS® Plus Instrument is an enhanced version of the CliniMACS Cell Selection Device. While the basic CliniMACS Instrument was designed to enrich stem and progenitor cells, the CliniMACS Plus Instrument provides an open platform technology enabling additional isolation strategies for many cell types, such as monocytes, dendritic cells, NK cells, T cells, and B cells (see chapter “CliniMACS Reagents” or call for application updates).

The combination of immunomagnetic reagent and tubing set offers high flexibility and versatility for individual separation strategies. Different-scale applications from rare-cell enrichment to bulk depletion are feasible.

Technical specifications

CliniMACS Plus Instrument

Model:	CS2 CE/UL
Size:	70 cm (w) × 60 cm (d), 90–130 cm (h)
Weight:	35 kg
Input voltage:	100–240 V AC
Power consumption:	180 VA

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.



CliniMACS® Plus Upgrade Kit “Software”

Product	Availability	Order no.
CliniMACS Plus Upgrade Kit “Software”	Europe, CAN, AUS, US ¹⁾	755-16
For availability in your country please contact your local representative.		
1) In the USA, the CliniMACS System components including the CliniMACS Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).		

Description

All CliniMACS® Plus Instruments can be upgraded to the latest status of the software.

The upgrade is CE-compliant.

Existing basic CliniMACS Instruments can be upgraded.

Kit components

CliniMACS Plus Upgrade Kit “Software” (755-16):

- Software Card
- CliniMACS User Manual
- 2 Bag Hanger assemblies
- Stabilization foot

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

Service contracts

CliniMACS® Service Contracts		
Product	Availability	Order no.
CliniMACS Full Service Agreement	Europe	160-000-973
CliniMACS Maintenance	Europe	160-000-974
CliniMACS Premium Service Agreement	CAN, US	160-000-701
CliniMACS Basic Service Agreement	CAN, US	160-000-702
For availability in your country please contact your local representative.		

CliniMACS® Tubing Sets

CliniMACS® Tubing Set and Tubing Set LS					
Product	Components	Regulatory status	Capacity (max.)	Availability	Order no.
CliniMACS Tubing Set	1 Tubing Set	CE	60×10 ⁹ total cells ²⁾	Europe, CAN, AUS, US ¹⁾	161-01
CliniMACS Tubing Set LS	1 Tubing Set	CE	120×10 ⁹ total cells ²⁾	Europe, CAN, AUS, US ¹⁾	162-01

For availability in your country please contact your local representative.

1) In the USA, the CliniMACS® System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

2) Capacity (max.) of labeled cells depends on the software program used.

Description

CliniMACS® Tubing Sets consist of pre-assembled tubing including a pre-column and a selection column. The fluid path of the tubing sets is sterile and non-pyrogenic. One package unit contains one sealed tubing set for single use.

The CliniMACS Tubing Set (161-01): The CliniMACS Tubing Set was designed for the separation of cells from up to 60×10⁹ total cells.

The CliniMACS Tubing Set LS (162-01): The CliniMACS Tubing Set LS was designed for the separation of cells from up to 120×10⁹ total cells.

Please inquire about recommended device and accessories.

Applications

The CliniMACS® Tubing Sets for clinical use were developed for the enrichment or depletion of cells from human heterogeneous hematologic cell populations in combination with the CliniMACS® System.

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

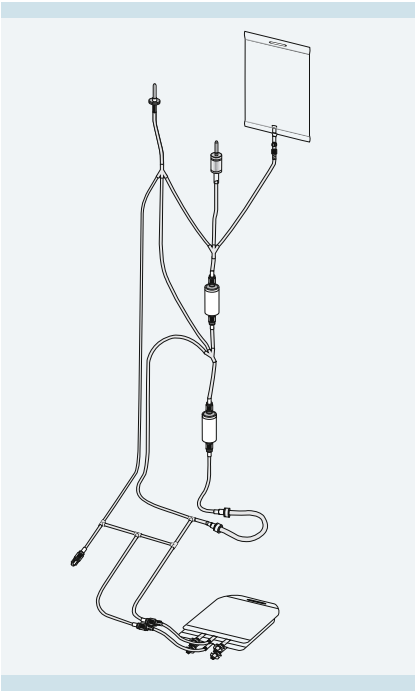


Figure 1: The CliniMACS Tubing Set.

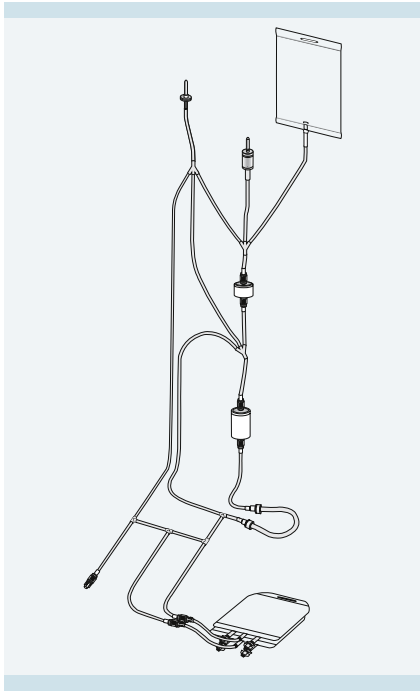


Figure 2: The CliniMACS Tubing Set LS.

CliniMACS® Tubing Sets for Research Use					
Product	Components	Regulatory status	Capacity (max.)	Availability	Order no.
CliniMACS Tubing Set for Research Use	1 Tubing Set	For research use only, not for use in humans ²⁾	60×10 ⁹ total cells ¹⁾	US	165-01
CliniMACS Tubing Set LS for Research Use	1 Tubing Set	For research use only, not for use in humans ²⁾	120×10 ⁹ total cells ¹⁾	US	168-01

For availability in your country please contact your local representative.
1) Capacity (max.) of labeled cells depends on the software program used.
2) Manufactured by Miltenyi Biotec GmbH under an ISO 9001 certified quality system.

Description

CliniMACS® Tubing Sets for Research Use consist of pre-assembled tubing including a pre-column and a selection column. The fluid path of the tubing sets is sterile and non-pyrogenic. One package unit contains one sealed tubing set for single use.

CliniMACS Tubing Set for Research Use (165-01):
The CliniMACS Tubing Set was designed for the separation of cells from up to 60×10⁹ total cells.

CliniMACS Tubing Set LS for Research Use (168-01):
The CliniMACS Tubing Set LS was designed for the separation of cells from up to 120×10⁹ total cells.

Please inquire about recommended device and accessories.

Applications

The CliniMACS® Tubing Sets for Research Use were developed for the enrichment or depletion of cells from human heterogeneous hematologic cell populations in combination with the CliniMACS® System.

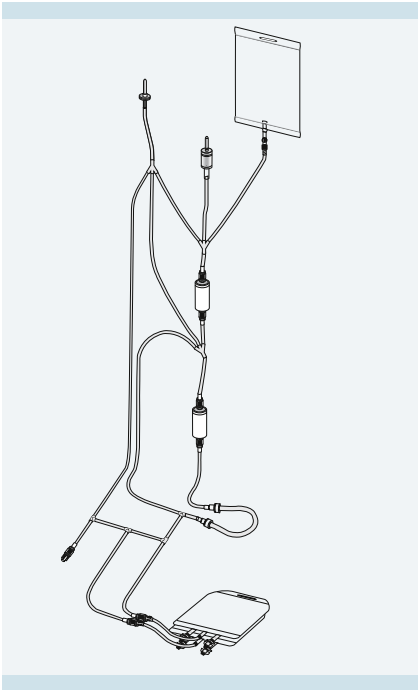


Figure 1: Schematic illustration of the CliniMACS Tubing Set for Research Use containing a pre-column and a selection column.

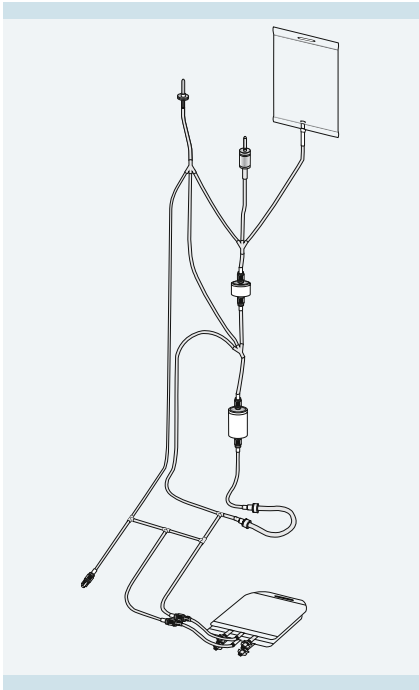


Figure 2: Schematic illustration of the CliniMACS Tubing Set LS for Research Use containing a pre-column and a selection column.

CliniMACS® Depletion Tubing Sets

Product	Components	Regulatory status	Capacity (max.)	Availability	Order no.
CliniMACS Depletion Tubing Set	1 Tubing Set	CE	40×10 ⁹ labeled cells from 120×10 ⁹ total cells	Europe, US ¹⁾	261-01
CliniMACS Depletion Tubing Set for Research Use	1 Tubing Set	For research use only, not for use in humans ²⁾	40×10 ⁹ labeled cells from 120×10 ⁹ total cells	US, CAN, AUS	266-01

For availability in your country please contact your local representative.

1) In the USA, the CliniMACS® System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

2) Manufactured by Miltenyi Biotec GmbH under an ISO 9001 certified quality system.

Description

CliniMACS® Depletion Tubing Sets contain tubing with a pre-assembled Cell Collection Bag, Reapplication Bag, Non-Target Cell Bag, and Buffer Waste Bag. The fluid path of the tubing sets is sterile and non-pyrogenic. One package unit contains one sealed tubing set for single use.

CliniMACS® Depletion Tubing Set (261-01):

The CliniMACS Depletion Tubing Set was designed for the depletion of up to 40×10⁹ labeled cells from up to 120×10⁹ total cells.

CliniMACS® Depletion Tubing Set for Research Use (266-01):

The CliniMACS Depletion Tubing Set for Research Use was designed for the depletion of up to 40×10⁹ labeled cells from up to 120×10⁹ total cells.

Please inquire about recommended device and accessories. The CliniMACS Plus Instrument with Software Version 2.40 or higher is required.

Applications

The CliniMACS Depletion Tubing Set is specifically designed for depletion of large quantities of cells from human heterogeneous hematologic cell populations in combination with the CliniMACS® System. It is also designed for a faster depletion of more cells as compared to other CliniMACS Tubing Sets.^{1,2}

References

1. Fritsch *et al.* (2004) Bone Marrow Transplant. 33 (Suppl.1): 595.
2. Fritsch *et al.* (2004) Cytotherapy 6(4): 432.

Related CliniMACS® Products

CliniMACS Plus Upgrade Kit "Software"	755-16
---------------------------------------	--------

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

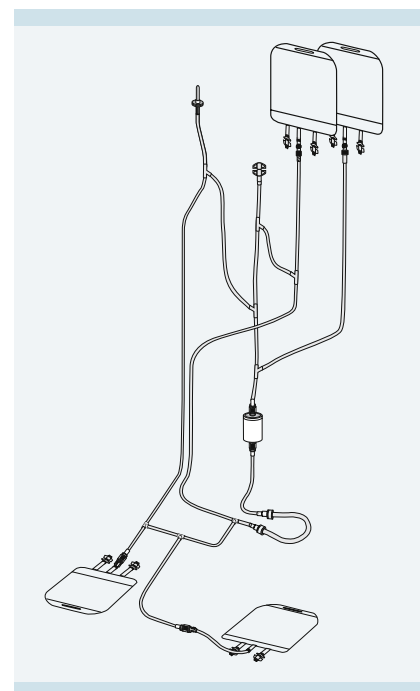


Figure 1: CliniMACS Depletion Tubing Set

CliniMACS® Tubing Sets 150 and 600 for Research Use					
Product	Components	Regulatory status	Capacity (max.)	Availability	Order no.
CliniMACS Tubing Set 600 for Research Use	5 Tubing Sets	For research use only, not for use in humans ¹⁾	20×10 ⁹ total cells	Worldwide	166-01
CliniMACS Tubing Set 150 for Research Use	5 Tubing Sets	For research use only, not for use in humans ¹⁾	10×10 ⁹ total cells	Worldwide	167-01

For availability in your country please contact your local representative.

1) Manufactured by Miltenyi Biotec GmbH under an ISO 9001 certified quality system.

Description

CliniMACS® Tubing Sets 600 and 150 for Research Use are ideal tools for pre-clinical cell separations on the CliniMACS Platform. The CliniMACS Tubing Sets for Research Use with reduced tubing geometry consist of a selection column without a pre-column. One package unit contains five separately sealed tubing sets for single use.

CliniMACS Tubing Set 600 for Research Use (166-01):

The CliniMACS Tubing Set 600 for Research Use was designed for the separation from up to 20×10⁹ total cells.

CliniMACS Tubing Set 150 for Research Use (167-01):

The CliniMACS Tubing Set 150 for Research Use was designed for the separation from up to 10×10⁹ total cells.

Please inquire about recommended device and accessories.

Applications

The CliniMACS Tubing Sets 600 and 150 for Research Use can be utilized for various pre-clinical separations. The system allows for large-scale, sterile cell separation. The cells can be used, for example, for animal models, cell culture assays, or other research-related applications.

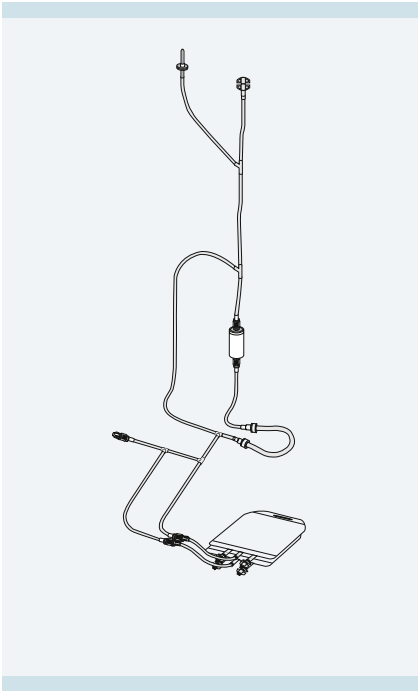


Figure 1: CliniMACS Tubing Set 600 for Research Use.

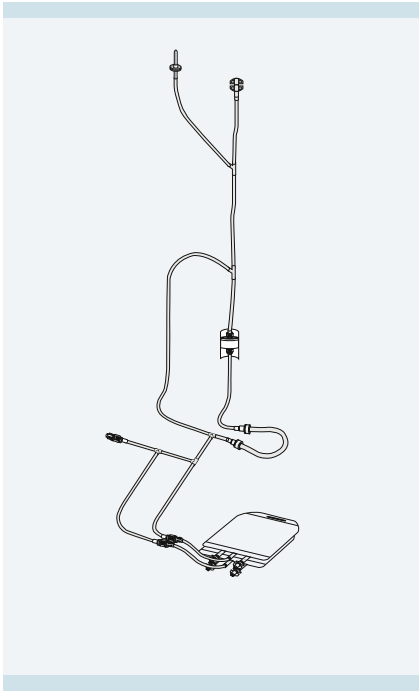


Figure 2: CliniMACS Tubing Set 150 for Research Use.

Stem and progenitor cells

CliniMACS® CD34 Enrichment

Background

The CD34 antigen is expressed on hematopoietic progenitor cells constituting a small subpopulation of bone marrow cells and peripheral blood cells. After severe damage, for example, after myeloablative conditioning, the hematopoietic system can be reconstituted by transplantation of allogeneic or autologous CD34⁺ hematopoietic progenitor cells. The concomitant removal of T cells and B cells is important especially for allogeneic grafts. CD34 enrichment of progenitor cells in the autologous setting may be used for removing contaminating tumor cells from the graft (purging).

Performance

Since its introduction, CliniMACS® CD34 System has been used in numerous clinical trials to provide CD34⁺ progenitor cells with high purity and yield in conjunction with excellent passive depletion of unwanted cells. A review of the performance in the allogeneic setting is shown in table 1.

Clinical results

The concomitant T and B cell depletion may reduce the incidence of graft versus host disease (GVHD) and EBV-induced lymphoproliferative disease in allogeneic transplantation settings. This has been suggested in the HLA identical setting with related^{1,2} as well as unrelated^{3,4} donors. It is thought to be crucial in the

haploidentical setting where, due to the high degree of HLA mismatch, GVHD is the major cause of transplantation-related mortality^{5,6}. It has also been shown that CliniMACS CD34 Enrichment is a means to provide tumor cell-depleted stem cell grafts in lymphoma and neuroblastoma patients^{7,8}. An emerging field for CD34-enriched cells has recently been seen in the depletion of autoreactive cells from stem cell grafts for diseases such as rheumatoid arthritis, systemic lupus erythematosus, multiple sclerosis, and systemic sclerosis⁹⁻¹².

References

1. Elmaagachi *et al.* (2003) *Blood* 101(2): 446–453.
2. Urbano-Ispizua *et al.* (2001) *Bone Marrow Transplant*. 28: 349–354.
3. Bornhaeuser *et al.* (2000) *Haematologica* 85: 839–847.
4. Lang *et al.* (2003) *Blood* 101: 1630–1636.
5. Handgretinger *et al.* (2003) *Pediatr Transplant*. 7 (Suppl. 3): 51–55.
6. Aversa *et al.* (2004) *Springer Semin Immunopathol*. 26 (1–2): 155–68.
7. Flohr *et al.* (2002) *Bone Marrow Transplant*. 29: 769–775.
8. Handgretinger *et al.* (2002) *Bone Marrow Transplant*. 29: 731–736.
9. Snowden *et al.* (2004) *J. Rheumatol* 31: 482–488.
10. Jayne *et al.* (2004) *Lupus* 13: 168–176.
11. Carreras *et al.* (2003) *Haematologica* 88: 306–314.
12. Rosen *et al.* (2000) *Arthritis Research* 2: 327–336.
13. Handgretinger *et al.* (2003) *Pediatr Transplant*. 7 (Suppl. 3): 51–55.
14. Lang *et al.* (2004) *ISCT Dublin, Miltenyi Biotec Satellite Symposium*.
15. Falzetti *et al.* (2002) *Bone Marrow Transplant*. 29 (Suppl. 2): S199.
16. Martin-Henao *et al.* (2002) *ISCT Barcelona*, #100.

	n = 32	n > 500	n = 293	n = 136	n = 30
Purity CD34 (%)	97.2	95–99	97.5	90	96
Recovery CD34 (%)	60	70	77	81	64
CD3 log depletion	4.6	> 5	4.6	4.6	5
CD19 log depletion	nd	< 0.5 %	nd	3.2	3.2

Table 1: CliniMACS CD34 Enrichment results from PBSCs of allogeneic donors (nd = not determined). References 1, 13–16.

CliniMACS® CD34 Cell Enrichment

Product	Regulatory status	Capacity (max.)	Availability	Order no.
CliniMACS CD34 Reagent	CE	0.6 × 10 ⁹ CD34 ⁺ cells from 60 × 10 ⁹ total cells	Europe, CAN, AUS, US ¹⁾	171-01
CliniMACS CD34 Complete Kit	CE-marked components	0.6 × 10 ⁹ CD34 ⁺ cells from 60 × 10 ⁹ total cells	Europe	178-01
CliniMACS CD34 Complete Kit with Accessories	CE-marked components	0.6 × 10 ⁹ CD34 ⁺ cells from 60 × 10 ⁹ total cells	Europe	177-01
CliniMACS CD34 Complete LS Kit	CE-marked components	1.2 × 10 ⁹ CD34 ⁺ cells from 120 × 10 ⁹ total cells	Europe	170-01
CliniMACS CD34 Complete LS Kit with Accessories	CE-marked components	1.2 × 10 ⁹ CD34 ⁺ cells from 120 × 10 ⁹ total cells	Europe	169-01
CliniMACS CD34 MicroBeads	For research use only, not for use in humans	0.6 × 10 ⁹ CD34 ⁺ cells from 60 × 10 ⁹ total cells	US	175-01

For availability in your country please contact your local representative.

1) In the USA, the CliniMACS® System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

Description

The CD34 antigen is a single transmembrane glycoprotein that is mainly expressed on human hematopoietic stem and progenitor

cells but is also present on endothelial progenitor cells. The CD34 antigen is involved in cell adhesion and is thought to function as a signaling molecule.

The CliniMACS® CD34 Products consist of superparamagnetic iron dextran particles directly conjugated to CD34 antibodies.

Stem and progenitor cells

One vial contains 7.5 mL sterile, non-pyrogenic solution suitable for magnetic labeling of up to 0.6×10^9 CD34⁺ cells from up to 60×10^9 total cells.

Please inquire about recommended device and accessories.

Applications

CliniMACS CD34 Products were developed for the enrichment of CD34⁺ cells from human heterogeneous hematologic cell populations in combination with the CliniMACS® System.


Referenced literature

CD34 enrichment of progenitor cells is performed to provide an allogeneic stem cell graft highly purified for CD34⁺ cells and depleted of unwanted cells.¹⁻¹¹ In autologous transplantation CD34⁺ progenitor cells can be enriched in order to passively remove tumor cells from the graft (purging).¹²⁻¹³

References

1. Elmaagacli *et al.* (2003) *Blood* 101(2): 446–453.
2. Bornhaeuser *et al.* (2000) *Haematologica* 85: 839–847.
3. Lang *et al.* (2003) *Blood* 101: 1630–1636.
4. Handgretinger *et al.* (2001) *Bone Marrow Transplant*. 27: 777–783.
5. Aversa *et al.* (2002) *Int. J. Hematol.* 76(Suppl.1): 165–168.
6. Chakraverty *et al.* (2002) *Blood* 99(3): 1071–1078.
7. Craddock *et al.* (2000) *Br. J. Haematol* 111(3): 797–800.
8. Gaipa *et al.* (2003) *Bone Marrow Transplant* 31: 857–860.
9. Lang *et al.* (2004) *Bone Marrow Transplant* 33: 25–32.
10. Benesch *et al.* (2004) *Br. J. Haematol* 125(1): 58–63.
11. Handgretinger *et al.* (2003) *Pediatr. Transplant*. 7(Suppl.3): 51–55.
12. Flohr *et al.* (2002) *Bone Marrow Transplant* 29: 769–775.
13. Handgretinger *et al.* (2002) *Bone Marrow Transplant* 29: 731–736.

A) Before enrichment

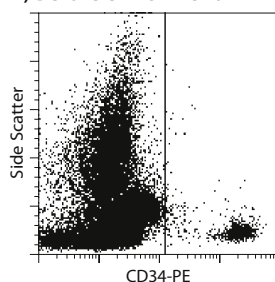

B) After enrichment

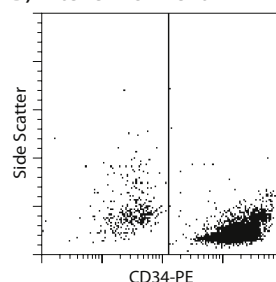


Figure 1: Flow cytometric analysis of CD34-enriched cells from mobilized leukapheresis product using the CliniMACS CD34 System. Cells are gated on viable CD45⁺ cells.

Kit components (not available in the US)

CliniMACS CD34 Complete Kit (178-01):

- 1 vial CliniMACS CD34 Reagent
- 1 CliniMACS Tubing Set
- 3 × 1 L CliniMACS PBS/EDTA Buffer

CliniMACS CD34 Complete Kit with Accessories (177-01):

- 1 vial CliniMACS CD34 Reagent
- 1 CliniMACS Tubing Set
- 3 × 1 L CliniMACS PBS/EDTA Buffer
- 1 Pre-System Filter
- 1 Luer/Spike Interconnector
- 1 Transfer Bag 150 mL

CliniMACS CD34 Complete LS Kit (170-01):

- 2 vials CliniMACS CD34 Reagent
- 1 CliniMACS Tubing Set LS
- 3 × 1 L CliniMACS PBS/EDTA Buffer

CliniMACS CD34 Complete LS Kit with Accessories (169-01):

- 2 vials CliniMACS CD34 Reagent
- 1 CliniMACS Tubing Set LS
- 3 × 1 L CliniMACS PBS/EDTA Buffer
- 1 Pre-System Filter
- 1 Luer/Spike Interconnector
- 1 Transfer Bag 150 mL

Related MACS® Products

CD34-FITC, human	130-081-001
CD34-PE, human	130-081-002
CD34-APC, human	130-090-954
CD133/2 (293C3)-PE, human	130-090-853
CD133/2 (293C3)-APC, human	130-090-854
CD45-FITC, human	130-080-202

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research only and not for therapeutic or diagnostic use.

Stem and progenitor cells

CliniMACS® CD133 (AC133) Enrichment

Background

The CD133 antigen has been identified as a stem cell marker. CD133 is specifically expressed on an immature subset of CD34⁺ cells, on circulating endothelial progenitor cells, as well as on a CD34⁺ stem cell subset (see figure 1)¹. CD133 is a marker for very early hematopoietic stem cells with high proliferative activity². The high potential for hematopoietic engraftment of isolated CD133⁺ cells has been shown in NOD/SCID repopulation assays^{3,4}. CD133-enriched cells have also been used as a starting fraction for the *ex vivo* expansion of hematopoietic progenitor cells from cord blood⁵⁻⁸. Clinically, CD133-enriched stem cell grafts have been employed in autologous^{9,10} and allogeneic transplantation, not only in the haploidentical^{11,12}, but also in the HLA-matched¹³ setting. Additionally, CD133⁺ stem cells can be utilized in non-hematological

applications such as regenerative medicine. CD133 is expressed on stem/progenitor cells of different tissues including bone marrow, peripheral blood, cord blood, and liver. CD133⁺ stem cells have been shown to harbor the capability to differentiate *in vitro* into cell types of various tissues.

Bone marrow-derived CD133⁺ stem cells have come into focus in non-hematological applications, especially ischemic heart diseases. Clinical trials are currently being conducted to address safety and feasibility of this approach¹⁴⁻¹⁸.

Performance

Since its introduction, the CliniMACS® CD133 System has demonstrated excellent performance providing CD133⁺ progenitor cells with high purity and yield in conjunction with excellent passive depletion of unwanted cells (see table 1).

References

1. De Wynter *et al.* (1998) *Stem Cells* 16: 387–396.
2. Goussetis *et al.* (2000) *J. Hematother. & Stem Cell Res.* 9: 827–840.
3. Kuci *et al.* (2003) *Blood* 101: 869–876.
4. Handgretinger *et al.* (2003) *Ann. N. Y. Acad. Sci.* 996: 141–151.
5. Pasino *et al.* (2000) *Brit. J. Haematol.* 108: 793–800.
6. Forraz *et al.* (2002) *Brit. J. Haematol.* 119: 516–524.
7. Kobari *et al.* (2001) *J. Hematother. & Stem Cell Res.* 10: 273–281.
8. Peled *et al.* (2004) *Cytother.* 6: 344–355.
9. Köhl *et al.* (2002) *Bone Marrow Transpl.* 29: 927–930.
10. Falzetti *et al.* (2002) *Blood* 100(11): 831a.
11. Lang *et al.* (2004) *Brit. J. Haematol.* 124: 72–79.
12. Lang *et al.* (2004) *Bone Marrow Transpl.* 33: 879–880.
13. Bornhaeuser *et al.* (2005) *Leukemia* 19: 161–165.
14. Pompilio *et al.* (2004) *Ann. Thorac. Surg.* 78(5): 1808–1812.
15. Klein *et al.* (2004) *The Heart Surgery Forum.* Vol. 7, Issue 5.
16. Ghodsizad *et al.* (2004) *Cytotherapy.* 6(5): 523–526.
17. Stamm *et al.* (2004) *Thorac. Cardiovasc. Surg.* 52(3): 152–158.
18. Stamm *et al.* (2003) *Lancet.* 361(9351): 45–46.
19. Gordon *et al.* (2003) *Bone Marrow Transpl.* 31: 17–22.
20. Lang *et al.* (2004) *ISCT Dublin, Miltenyi Biotec Satellite Symposium.*
21. Bornhaeuser *et al.* (2005) *Leukemia* 19: 161–165.

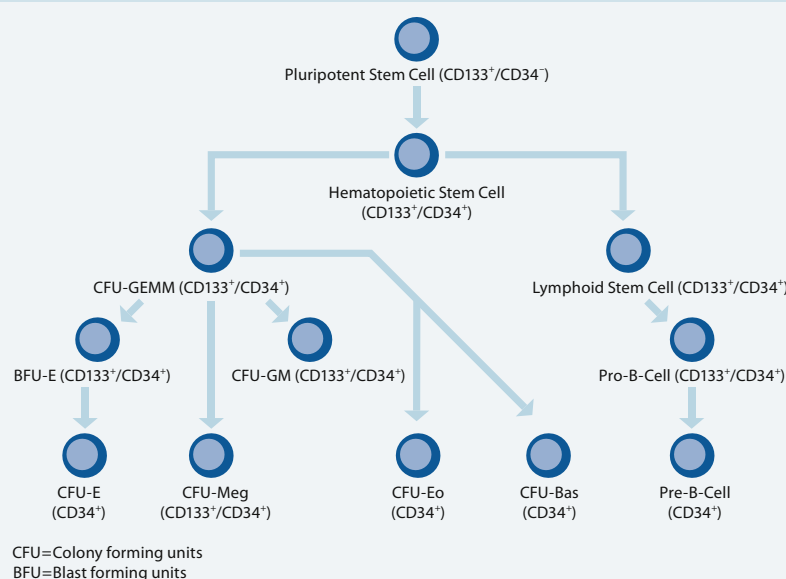


Figure 1: The early hematopoietic progenitor cell marker CD133 is ancestral to CD34.

	n = 11	n = 84	n = 12
Purity CD133 (%)	94	93	97
Recovery CD133 (%)	69	81	69
CD3 log depletion	4.2	3.8	nd

Table 1: CliniMACS CD133 enrichment results from PBSCs of allogeneic donors (nd=not determined). References 19–21.

CliniMACS® CD133 (AC133) Cell Enrichment

Product	Regulatory status	Capacity (max.)	Availability	Order no.
CliniMACS CD133 Reagent	CE	0.6×10^9 CD133 ⁺ cells from 60×10^9 total cells	Europe, US ¹⁾	172-01
CliniMACS CD133 Complete Kit	CE-marked components	0.6×10^9 CD133 ⁺ cells from 60×10^9 total cells	Europe	196-01
CliniMACS CD133 Complete Kit with Accessories	CE-marked components	0.6×10^9 CD133 ⁺ cells from 60×10^9 total cells	Europe	198-01
CliniMACS CD133 Complete LS Kit	CE-marked components	1.2×10^9 CD133 ⁺ cells from 120×10^9 total cells	Europe	197-01
CliniMACS CD133 Complete LS Kit with Accessories	CE-marked components	1.2×10^9 CD133 ⁺ cells from 120×10^9 total cells	Europe	199-01
CliniMACS CD133 MicroBeads	For research use only, not for use in humans	0.6×10^9 CD133 ⁺ cells from 60×10^9 total cells	US, CAN, AUS	195-01

For availability in your country please contact your local representative.

1) In the USA, the CliniMACS® System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

Description

CD133 is expressed on immature hematopoietic stem and progenitor cells but is not found on mature blood cells. In contrast to the CD34 antigen, CD133 is not expressed by late progenitors such as pre-B cells, CFU-E, and CFU-G. The stem cell marker CD133 was formerly known as AC133.

CliniMACS® CD133 Products consist of superparamagnetic iron dextran particles directly conjugated to CD133 antibodies.

One vial contains 7.5 mL sterile, non-pyrogenic solution suitable for magnetic labeling of up to 0.6×10^9 CD133⁺ cells from up to 60×10^9 total cells. Please inquire about recommended device and accessories.


Applications

CliniMACS CD133 (AC133) Products were developed for the enrichment of CD133⁺ cells from human heterogeneous hematologic cell populations in combination with the CliniMACS® System.

Referenced literature

CD133 enrichment of progenitor cells is performed to provide an allogeneic stem cell graft highly purified for CD133⁺ cells and depleted of unwanted cells.¹⁻⁶ In autologous transplantation CD133⁺ progenitor cells can be enriched in order

A) Before enrichment

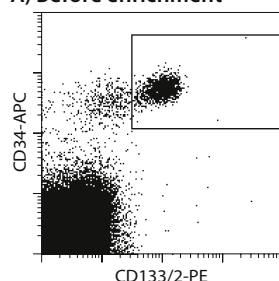

B) After enrichment


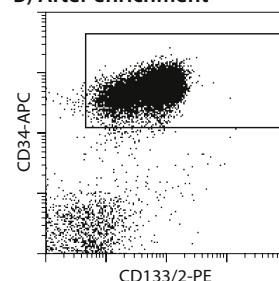


Figure 1: Flow cytometric analysis of CD133-enriched cells from a mobilized leukapheresis product using the CliniMACS CD133 System.

A) Before enrichment

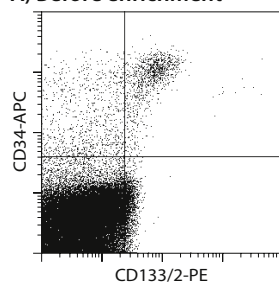

B) After enrichment

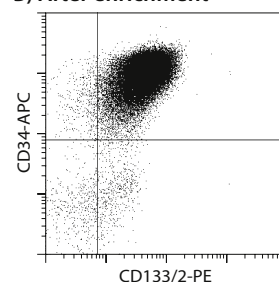


Figure 2: Flow cytometric analysis of CD133-enriched stem cells from a bone marrow aspirate using the CliniMACS CD133 System.

to passively remove tumor cells from the graft (purging).⁷ In addition, stem cells enriched for the CD133 antigen are being explored for use in regenerative medicine (e.g. cardiac diseases).⁸⁻¹²

References

1. Bitan *et al.* (2005) *Exp. Hematol.* 33: 713–718.
2. Bornhäuser *et al.* (2005) *Leukemia* 19: 161–165.
3. Peled *et al.* (2004) *Cytotherapy* 6: 344–355.
4. Lang *et al.* (2004) *Brit. J. Haematol.* 124: 72–79.
5. Lang *et al.* (2004) *Bone Marrow Transplant.* 33 (1): 879–880.
6. Gordon *et al.* (2003) *Bone Marrow Transplant.* 31: 17–22.
7. Koehl *et al.* (2002) *Bone Marrow Transplant.* 29: 927–930.
8. Pompilio *et al.* (2004) *Ann. Thorac. Surg.* 78 (5): 1808–1812.
9. Klein *et al.* (2004) *The Heart Surgery Forum.* Vol. 7, Issue 5.
10. Ghodsizad *et al.* (2004) *Cytotherapy* 6 (5): 523–526.
11. Stamm *et al.* (2004) *Thorac Cardiovasc. Surg.* 52 (3): 152–158.
12. Stamm *et al.* (2003) *Lancet* 361 (9351): 45–46.

Stem and progenitor cells

Kit components (not available in the US)

CliniMACS CD133 Complete Kit

(196-01):

- 1 vial CliniMACS CD133 Reagent
- 1 CliniMACS Tubing Set
- 3×1 L CliniMACS PBS/EDTA Buffer

CliniMACS CD133 Complete Kit with

Accessories (198-01):

- 1 vial CliniMACS CD133 Reagent
- 1 CliniMACS Tubing Set
- 3×1 L CliniMACS PBS/EDTA Buffer
- 1 Pre-System Filter
- 1 Luer/Spike Interconnector
- 1 Transfer Bag 150 mL

CliniMACS CD133 Complete LS Kit

(197-01):

- 2 vials CliniMACS CD133 Reagent
- 1 CliniMACS Tubing Set LS
- 3×1 L CliniMACS PBS/EDTA Buffer

CliniMACS CD133 Complete LS Kit with

Accessories (199-01):

- 2 vials CliniMACS CD133 Reagent
- 1 CliniMACS Tubing Set LS
- 3×1 L CliniMACS PBS/EDTA Buffer
- 1 Pre-System Filter
- 1 Luer/Spike Interconnector
- 1 Transfer Bag 150 mL

Related MACS® Products

CD133/2 (293C3)-PE, human	130-090-853
CD133/2 (293C3)-APC, human	130-090-854
CD34-FITC, human	130-081-001
CD34-PE, human	130-081-002
CD34-APC, human	130-090-954
CD45-FITC, human	130-080-202
CD45-PE, human	130-080-201
CD45-APC, human	130-091-230

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research only and not for therapeutic or diagnostic use.

CliniMACS® T and B cell depletion (CD3/CD19 combination)¹⁾

Using the CliniMACS® System together with the CliniMACS CD3/CD19 Combination makes it possible to simultaneously deplete T cells and B cells from a graft.

In contrast to stem cell enrichment, this method preserves NK cells and other effector cells in the graft (see also table 1).¹

Safety and feasibility of CD3/CD19-depleted grafts in the haploidentical and unrelated donor setting have been explored in a number of clinical studies²⁻⁵.

Table 1 summarizes the CD3/CD19 depletion results with the CliniMACS System and the CliniMACS Tubing Set LS.

However, with the CliniMACS Tubing Set LS the depletion procedure is limited to a maximum of 20×10^9 labeled cells. Therefore, a new tubing set was developed, called the CliniMACS Depletion Tubing Set, which has been designed for depletion of up to 40×10^9 labeled cells with a reduced processing time. The new procedure uses the program DEPLETION 3.1 (software version 2.40 or higher).

	n = 5	n = 25	n = 6	n = 14
Cell source	PBSCs	PBSCs	PBSCs	BM
Recovery CD34 (%)	69	70	77	67
Recovery NK cells (%)	36	nd	67	75
Remaining T cells (%)	0.02	0.01	nd	0.11
Remaining B cells (%)	0.1	0.005	nd	1.6
CD3 log depletion	3.4	3.9	3.7	3
CD19 log depletion	2.2	nd	3.7	nd

Table 1: CliniMACS CD3/CD19 Depletion results from allogeneic settings (nd=not determined). References 1, 3–5.

References

1. Barfield *et al.* (2004) *Cytother* 6: 1–6.
2. Benaim *et al.* (2003) *Blood* 102(11): 969a.
3. Lang *et al.* (2004) ISCT Dublin, Miltenyi Biotec Satellite Symposium.
4. Fritsch *et al.* (2004) ISCT Dublin, Ireland #192.
5. Preijers *et al.* (2004) *Bone Marrow Transplant*. 33 (Suppl 1): S93.

1) Not available in the United States.

CliniMACS® CD3/CD19 Depletion

Product	Regulatory status	Capacity (max.)	Availability	Order no.
CliniMACS CD3/CD19 Reagent Combination	CE-marked components	5×10^9 CD19 ⁺ cells and 15×10^9 CD3 ⁺ cells from 40×10^9 total cells	Europe	704-00
CliniMACS CD3/CD19 Complete Kit	CE-marked components	5×10^9 CD19 ⁺ cells and 15×10^9 CD3 ⁺ cells from 40×10^9 total cells	Europe	740-01
CliniMACS CD3/CD19 Large-Scale Complete Kit	CE-marked components	10×10^9 CD19 ⁺ cells and 30×10^9 CD3 ⁺ cells from 80×10^9 total cells	Europe	740-02

For availability in your country please contact your local representative

Description

The CliniMACS® CD3/CD19 Products consist of superparamagnetic iron dextran particles directly conjugated to CD3 and CD19 antibodies.

Each vial contains 7.5 mL sterile, non-pyrogenic solution suitable for magnetic labeling of up to 15×10^9 CD3⁺ cells and 5×10^9 CD19⁺ cells from up to 40×10^9 total cells, respectively.

Please inquire about recommended device and accessories.

Applications

The CliniMACS® CD3/CD19 Products were developed for the simultaneous depletion of unwanted T and B cells from human heterogeneous hematologic cell populations in combination with the CliniMACS® System. This approach keeps stem and progenitor cells untouched and leaves immune effector cells, such as NK cells and dendritic cells, in the cellular product.¹⁻²

Kit components (not available in the US)

CliniMACS CD3/CD19 Reagent Combination (704-00):

- 1 vial CliniMACS CD3 Reagent
- 1 vial CliniMACS CD19 Reagent


CliniMACS CD3/CD19 Complete Kit (740-01):

- 1 vial CliniMACS CD3 Reagent
- 1 vial CliniMACS CD19 Reagent
- 1 CliniMACS Depletion Tubing Set
- 3 × 1 L CliniMACS PBS/EDTA Buffer


CliniMACS CD3/CD19 Large-Scale Complete Kit (740-02):

- 2 vials CliniMACS CD3 Reagent
- 2 vials CliniMACS CD19 Reagent
- 1 CliniMACS Depletion Tubing Set
- 3 × 1 L CliniMACS PBS/EDTA Buffer

A) Before depletion

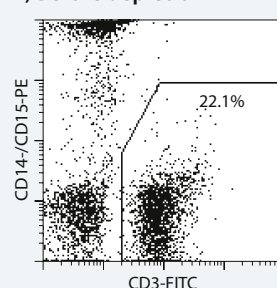


B) After depletion

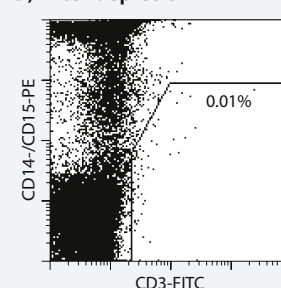


T cell depletion

C) Before depletion

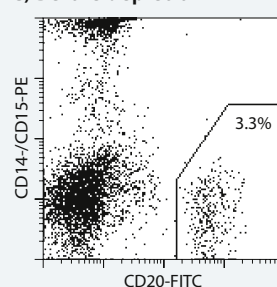

D) After depletion

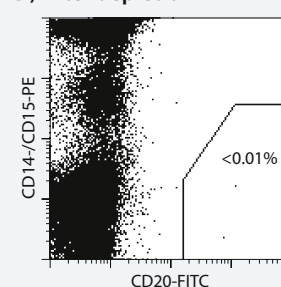

B cell depletion

Figure 1: Flow cytometric analysis of simultaneous depletion of T and B cells from a buffy coat using the CliniMACS CD3/CD19 System.

Related CliniMACS® Products

CliniMACS CD3/CD56 Reagent Combination ¹⁾	704-01
CliniMACS CD3 Reagent	273-01
CliniMACS CD19 Reagent	179-01
CliniMACS CD56 Reagent	271-01

1) Not available in the US.

Related MACS® Products

CD14-PE, human	130-091-242
CD15-PE, human	130-091-375
CD20-FITC, human	130-091-108
CD45-FITC, human	130-080-202

References

1. Barfield *et al.* (2004) *Cytotherapy* 6: 1–6.
2. Gordon *et al.* (2002) *Bone Marrow Transplant*. 30: 69–74.

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research only and not for therapeutic or diagnostic use.

CliniMACS® T Cell Separation

T cells play critical roles in the regulation of immune responses and are responsible for mediating both beneficial and harmful immune responses.

In the transplant setting, graft versus leukemia (GVL) effects and graft versus host diseases (GVHD) are in a large part conveyed by donor T lymphocytes as part of the unfolding adaptive immune response. New immunotherapy strategies have evolved, designed to separate GVL from GVHD, as well as to boost GVL and reduce GVHD.

The CliniMACS® System, together with the broadened CliniMACS T Cell Product portfolio, allows for enrichment or depletion of different T cell subsets providing the basis for a variety of different applications.

The CliniMACS CD3 Reagent was developed for the depletion of T cells from unmobilized as well as mobilized leukapheresis products. The rationale for this application is that the depletion of T cells may lower the risk of GVHD. Combined depletion of T and B cells from

the graft using CliniMACS CD3 and CD19 Reagents leaves stem and progenitor cells and, in addition, provides a graft which contains NK cells¹. Enrichment of NK cells can also be achieved using the CliniMACS CD56 Reagent after prior depletion of T cells with the CliniMACS CD3 Reagent.² For certain oncological settings preliminary data suggest that T cells in a graft or as a donor lymphocyte infusion (DLI) may be beneficial.³ In these settings a CD4 enrichment or CD8 depletion may be desired. CD4⁺ T cells not only specifically help CD8⁺ T cells during an immune reaction but also seem to support their prolonged survival in the recipient⁴. The depletion of CD8⁺ T cells from DLI is described in various protocols for the prevention of GVHD.⁵ The CliniMACS CD4 Reagent and CD8 Reagent are the reagents of choice for investigating potential applications based on such rationale. CD25 was first described as a T cell proliferation marker. It is now regarded as a marker which defines activated/reactive T cells.

Alloreactive T cells can cause severe complications (GVHD) in transplantation.⁶ Thus, it may be desirable to deplete CD25⁺ T cells from apheresis products. CD25^{hi}CD4⁺ regulatory T cells are the focus of T cell research as they have been shown to be capable of specifically modulating immune reactions. For a variety of applications, these modulating capacities could potentially be utilized.⁷ The CD25⁺ T cells can be depleted or enriched using the CliniMACS CD25 System.^{8,9}

Antigen-specific T cells have become another focus in T cell research and opportunities for utilizing such cells for clinical-scale applications are being exploited. In order for an antigen-specific CD8⁺ cell to survive long enough and thus be able to fight an infection or a tumor, the help of CD4⁺ T cells is mandatory. The CliniMACS Cytokine Capture System (IFN-gamma) is the first product developed that allows the concomitant clinical-scale enrichment of both CD4⁺ and CD8⁺ T cells specific for a given antigen.¹⁰

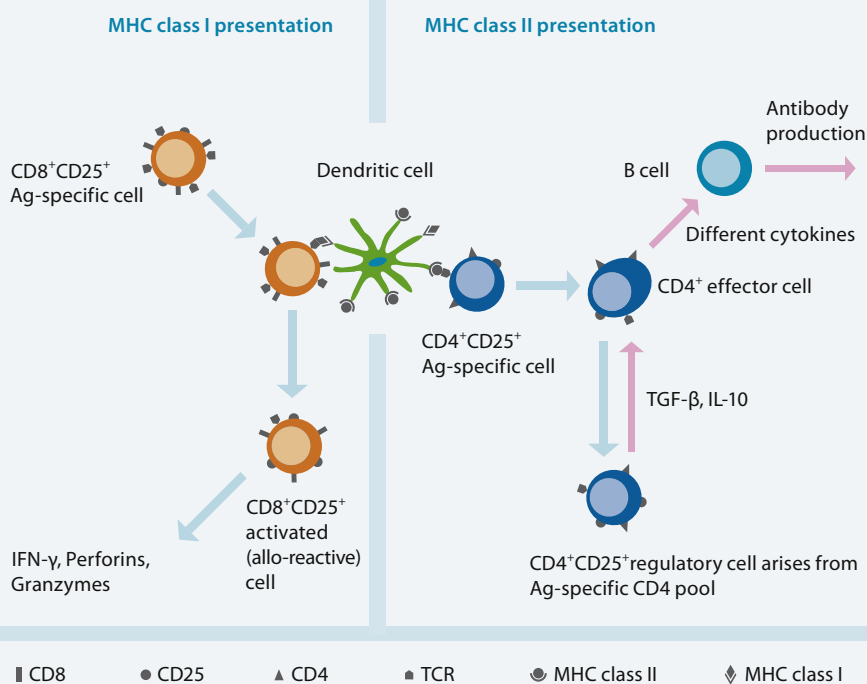


Figure 1: Schematic diagram of the interface between dendritic cells and T cells.

References

1. Barfield *et al.* (2004) *Cytotherapy* 6: 1–6.
2. Iyengar *et al.* (2003) *Cytotherapy* 5 (6): 479–484.
3. Falkenburg *et al.* (1999) *Blood* 94 (4): 1201–1208.
4. Rosenberg *et al.* (2004) *PNAS, USA* 101: 14639–14645.
5. Soiffer *et al.* (2002) *Biol. Blood Marrow Transplant* 8: 625–632.
6. André-Schmutz *et al.* (2002) *Lancet* 360: 130–137.
7. Hoffmann *et al.* (2004) *Blood* 104: 895–903.
8. Hoffmann *et al.* (2004) *Blood* 104: 2851.
9. Powell *et al.* (2005) *J. Immunother.* 28: 403–411.
10. Feuchtinger *et al.* (2004) *Exp. Hem.* 32: 282–289.

CliniMACS® CD3 Cell Depletion

Product	Regulatory status	Capacity (max.)	Availability	Order no.
CliniMACS CD3 Reagent	CE	15 × 10 ⁹ CD3 ⁺ cells from 40 × 10 ⁹ total cell	Europe, US ¹⁾	273-01
CliniMACS CD3 MicroBeads	For research use only, not for use in humans	15 × 10 ⁹ CD3 ⁺ cells from 40 × 10 ⁹ total cells	US, CAN, AUS	176-01

For availability in your country please contact your local representative.

1) In the USA, the CliniMACS® System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

Description

CD3 is expressed on all T cells and is associated with the T cell receptor. Approximately 70–80% of human peripheral blood lymphocytes and 65–85% of thymocytes are CD3⁺.

CliniMACS® CD3 Products consist of superparamagnetic iron dextran particles directly conjugated to CD3 antibodies.

One vial contains 7.5 mL sterile, non-pyrogenic solution suitable for magnetic labeling of up to 15 × 10⁹ CD3⁺ cells from up to 40 × 10⁹ total cells.

Please inquire about recommended device and accessories.


Applications

CliniMACS T Cell CD3 Products were developed for the depletion of unwanted T cells from human heterogeneous hematologic cell populations in combination with the CliniMACS® System. This approach maintains the stem cells and immune effector cells, such as NK cells, in the cellular product.^{1,2}

References

- Gordon *et al.* (2002) Bone Marrow Transplant. 30: 69–74
- Chen *et al.* (2006) Brit. J. Haem. 135: 524–532.

A) Before depletion

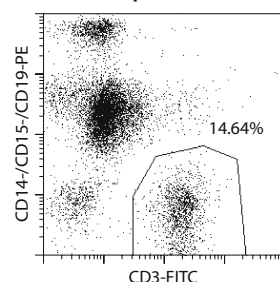

B) After depletion

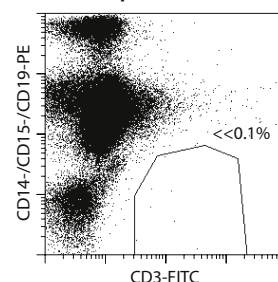


Figure 1: Flow cytometric analysis of CD3-depleted cells from a leukapheresis product using the CliniMACS CD3 System.

Related CliniMACS® Products

CliniMACS CD19 Reagent	179-01
CliniMACS CD56 Reagent	271-01
CliniMACS CD3/CD19 Reagent Combination ¹⁾	704-00

1) Not available in the US.

Related MACS® Products

CD14-PE, human	130-091-242
CD15-PE, human	130-091-375
CD19-PE, human	130-091-247

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research only and not for therapeutic or diagnostic use.

T cells

CliniMACS® CD4 Cell Separation

Product	Regulatory status	Capacity (max.)	Availability	Order no.
CliniMACS CD4 Reagent	CE	12×10 ⁹ CD4 ⁺ cells from 40×10 ⁹ total cells	Europe, US ¹⁾	276-01
CliniMACS CD4 MicroBeads	For research use only, not for use in humans	12×10 ⁹ CD4 ⁺ cells from 40×10 ⁹ total cells	US, CAN, AUS	304-01

For availability in your country please contact your local representative.

1) In the USA, the CliniMACS® System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

Description

CD4 is an accessory molecule in the recognition of foreign antigens in association with MHC class II antigens by T cells. T helper cells and to a lower degree monocytes and dendritic cells express CD4.

The CliniMACS® CD4 Products consist of superparamagnetic iron dextran particles directly conjugated to CD4 antibodies.

One vial contains 7.5 mL sterile, non-pyrogenic solution suitable for magnetic labeling of up to 12×10⁹ CD4⁺ cells from up to 40×10⁹ total cells.

Please inquire about recommended device and accessories.

Applications


CliniMACS® CD4 Products were developed for the depletion or enrichment of CD4⁺ cells from human heterogeneous hematologic cell populations in combination with the CliniMACS® System.

Referenced literature

In the allogeneic transplantation setting, CD4⁺ T cells in donor lymphocyte infusions may be capable of mounting anti-tumor responses, concomitantly reducing the risk of GVHD¹ and reconstituting the recipient's immune system.²

Recently, Th2 cells have been successfully used to support stem cell grafts.³ In addition, the use of CD4⁺ T cells may be the basis for various gene-therapeutic approaches such as in the HIV setting.⁴⁻⁶

A) Before enrichment

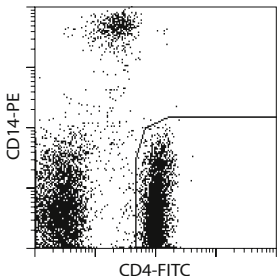

B) After enrichment


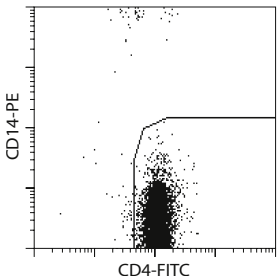


Figure 1: Flow cytometric analysis of CD4-enriched cells using the CliniMACS CD4 System.

A) Before depletion

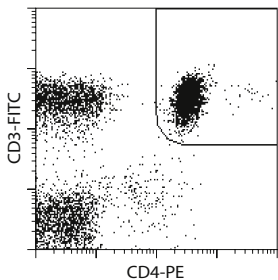

B) After depletion

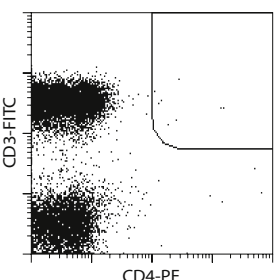


Figure 2: Flow cytometric analysis of CD4-depleted cells using the CliniMACS CD4 System.

References

1. Alyea *et al.* (1998) Blood 91: 3671–3680.
2. Bellucci *et al.* (2002) Blood 15: 4610–4617.
3. Fowler *et al.* (2004) Semin. Oncol. 31: 56–67.
4. Bernstein *et al.* (2004) Clin. Immunol. III: 262–274.
5. Levine *et al.* (2002) Nat. Med. 8: 47–53.
6. Davis *et al.* (2004) Mol. Ther. 9: 160–172.

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research only and not for therapeutic or diagnostic use.

Related MACS® Products

CD4-FITC, human	130-080-501
CD4-PE, human	130-091-231
CD4-APC, human	130-091-232
CD3-FITC, human	130-080-401
CD25-Biotin, human	130-091-235
Anti-Biotin-PE	130-090-756
Anti-Biotin-APC	130-090-856
CD8-FITC, human	130-080-601
CD30-PE, human	130-081-401
CD14-PE, human	130-091-242
CD25-PE, human	130-091-024
CD8-PE, human	130-091-084
CD8-APC, human	130-091-076
CD14-FITC, human	130-080-701
CD14-APC, human	130-091-243
Anti-Biotin-FITC	130-090-857

CliniMACS® CD8 Cell Depletion

Product	Regulatory status	Capacity (max.)	Availability	Order no.
CliniMACS CD8 Reagent	CE	4 × 10 ⁹ CD8 ⁺ cells from 40 × 10 ⁹ total cells	Europe, US ¹⁾	275-01
CliniMACS CD8 MicroBeads	For research use only, not for use in humans	4 × 10 ⁹ CD8 ⁺ cells from 40 × 10 ⁹ total cells	US, CAN, AUS	308-01

For availability in your country please contact your local representative.

1) In the USA, the CliniMACS® System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

Description

The CD8 antigen, a co-receptor for MHC class I molecules, is expressed strongly on cytotoxic T cells and dimly on a subset of NK cells.

The CliniMACS® CD8 Products consist of superparamagnetic iron dextran particles directly conjugated to CD8 antibodies.

One vial contains 7.5 mL sterile, non-pyrogenic solution suitable for magnetic labeling of up to 4 × 10⁹ CD8⁺ cells from up to 40 × 10⁹ total cells.

Please inquire about recommended device and accessories.

Applications

CliniMACS CD8 Products were developed for depletion of CD8⁺ T cells from human heterogeneous hematologic cell populations in combination with the CliniMACS System.


Referenced literature

In the allogeneic transplantation setting CD8⁺ T cells may be removed from a graft or DLI with the goal of preventing GVHD while maintaining GVL effects.¹⁻⁴ Additionally, the depletion of CD8⁺ T cells might be part of a strategy for the enrichment of regulatory T cells. Enrichment of CD8⁺ T cells is also an option for the generation of specific T cell lines.⁵

References

1. Nimer *et al.* (1994) *Transpl.* 57: 82–87.
2. Baron *et al.* (2003) *Haematologica* 88: 835–837.
3. Meyer *et al.* (2004) *Blood* 104: 5069.
4. Meyer *et al.* (2007) *Blood* 109: 374–382.
5. Savoldo *et al.* (2002) *Blood* 100: 4059–4066.

A) Before depletion

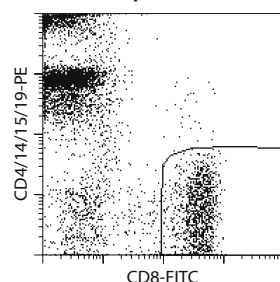

B) After depletion

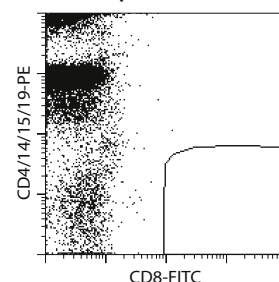


Figure 1: Flow cytometric analysis of CD8-depleted cells using the CliniMACS CD8 System.

Related MACS® Products

CD8-FITC, human	130-080-601
CD8-PE, human	130-091-084
CD8-APC, human	130-091-076
CD4-PE, human	130-091-231
CD4-APC, human	130-091-232
CD3-FITC, human	130-080-401
CD25-Biotin, human	130-091-235
Anti-Biotin-PE	130-090-756
Anti-Biotin-APC	130-090-856
CD30-PE, human	130-081-401
CD4-FITC, human	130-080-501

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research only and not for therapeutic or diagnostic use.

T cells

CliniMACS® CD25 Cell Separation

Product	Regulatory status	Capacity (max.)	Availability		Order no.
CliniMACS CD25 Reagent	CE	6 × 10 ⁹ CD25 ⁺ cells (highly expressing CD25) from 40 × 10 ⁹ total cells	Europe, US ¹⁾	NEW	274-01
CliniMACS CD25 MicroBeads	For research use only, not for use in humans	6 × 10 ⁹ CD25 ⁺ cells (highly expressing CD25) from 40 × 10 ⁹ total cells	US, CAN, AUS		325-01

For availability in your country please contact your local representative.

1) In the USA, the CliniMACS® System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

Description

The CD25 antigen, the low affinity interleukin-2 receptor alpha chain (IL-2R α), is expressed on activated T and B cells. In addition, CD25 is highly expressed on naturally occurring CD25^{hi}CD4⁺ regulatory T cells.

The CliniMACS® CD25 Products consist of superparamagnetic iron dextran particles directly conjugated to CD25 antibodies.

One vial contains 7.5 mL sterile, non-pyrogenic solution suitable for magnetic labeling of 6 × 10⁹ CD25⁺ cells (highly expressing CD25) from up to 40 × 10⁹ total cells.

Please inquire about recommended device and accessories.

Applications


CliniMACS® CD25 Products were developed for the enrichment or depletion of CD25⁺ cells from human heterogeneous hematologic cell populations in combination with the CliniMACS® System.

Referenced literature

In the allogeneic transplantation setting a promising approach to reduce risk of severe GVHD may be the selective removal of CD25⁺ alloreactive T cells.¹⁻³

Studies in the mouse model paved the way to investigate the role of CD25^{hi}CD4⁺ regulatory T cells for the improvement of stem cell transplantation. The addition of regulatory T cells, either to the graft or DLI may reduce the risk of GVHD.⁴ CD25^{hi}CD4⁺ regulatory T cells may also be

A) Before enrichment

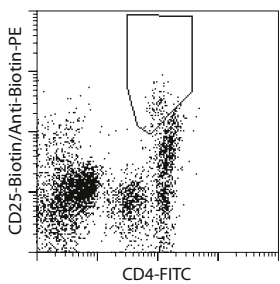

B) After enrichment


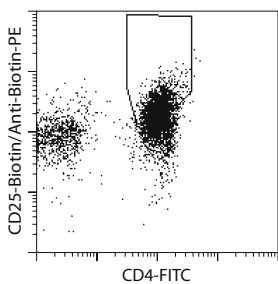


Figure 1: Flow cytometric analysis of CD25-enriched cells using the CliniMACS CD25 System.

Analysis of the content of CD25^{hi}CD4⁺ regulatory T cells. For flow cytometric analysis a specific electronic window is defined. Regulatory T cells have been described as CD4⁺ T cells which highly express CD25. The expression of CD4 is slightly reduced as compared to conventional CD4⁺ T helper cells.⁶

A) Before depletion

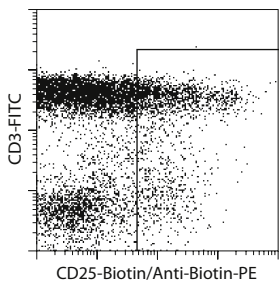

B) After depletion

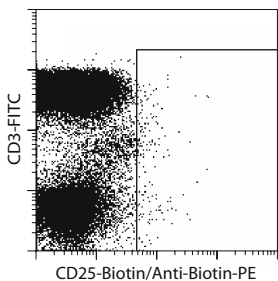


Figure 2: Flow cytometric analysis of CD25-depleted cells using the CliniMACS CD25 System.

depleted from patient's leukapheresis samples for immunotherapeutic approaches designed to favor anti-tumor responses.⁵

References

- Solomon *et al.* (2002) *Cytotherapy* 4 (5): 395–406.
- André-Schmutz *et al.* (2003) *Cytotherapy* 7 (2): 102–108.
- Amrolia *et al.* (2003) *Blood* 102 (6): 2292–2299.
- Edinger *et al.* (2003) *Nat. Med.* 9: 1144–1150.
- Powell *et al.* (2005) *J. Immunother.* 28(4): 403–11.
- Baecher-Allen *et al.* (2001) *J. Immunol.* 167: 1245–1253.

Related MACS® Products

CD25-Biotin, human	130-091-235
Anti-Biotin-PE	130-090-756
Anti-Biotin-APC	130-090-856
CD3-FITC, human	130-080-401
CD4-FITC, human	130-080-501
CD8-FITC, human	130-080-601

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research only and not for therapeutic or diagnostic use.

Antigen-specific T cells

CliniMACS® Antigen-specific T cell enrichment

Cytokine secretion is a hallmark of functionality of antigen-specific T cells. Large amounts of cytokines are produced by effector and memory T cells after short-term restimulation with “their” specific antigen.

For a multitude of challenges, it is of great interest to use antigen-specific T cells. Such applications could include prevention or managing of infection with Cytomegalovirus (CMV), Adenovirus, or Epstein-Barr Virus (EBV)^{1,2}. After immunosuppression (i.e. in the course of chemotherapy or organ-transplantation) such infections pose life-threatening complications, which sometimes are hard to fight by conventional means.

The treatment of tumors, which are resistant to conventional therapy, may become another field of application for antigen-specific T cells. Possible scenarios may also include autoimmune diseases such as rheumatoid arthritis and others.

It has been shown that optimal T cell immunity requires both CD4⁺ helper T cells and CD8⁺ cytotoxic T cells.³ The CliniMACS® Cytokine Capture System (IFN-gamma) enables the enrichment of viable antigen-specific CD4⁺ and antigen-specific CD8⁺ T cells based on the cytokines they secrete.^{4–6}

The CliniMACS Cytokine Capture System (IFN-gamma) offers unique reagents for clinical-scale enrichment of antigen-specific T cells, specifically developed for use exclusively on the CliniMACS® System.

Those cell preparations are the prerequisite for detailed analyses on the frequency and phenotype of the key players in a given immune response. Since the cells are viable after enrichment, subsequent *in vitro* steps such as expansion of the antigen-specific cells can be performed. Such an expansion step can also be performed “in the bag” to provide a greater degree of assurance of sterility.

The sensitivity and reliability of the assays have been demonstrated in a large panel of high-level scientific publications. These describe a variety of research applications from oncology or virology to autoimmune diseases.^{4–6}

Using the CliniMACS Cytokine Capture System (IFN-gamma) the enrichment can be performed directly “in the bag” starting from up to 1×10^9 total cells. As for all CliniMACS applications, the entire process of cell enrichment is performed in a closed, sterile system. The resulting viable cells may be directly used in subsequent steps of the respective protocol. Alternatively, the cells may be manipulated or expanded *in vitro* or phenotypically analyzed.

References:

1. Cohen *et al.* (2002) *Virology* 304: 474.
2. Bissinger *et al.* (2002) *Exp. Hematol.* 30: 1178.
3. Dudley *et al.* (2002) *Science* 298: 850.
4. Campbell (2003) *Methods* 31: 150–159.
5. Feuchtinger *et al.* (2004) *Exp. Hematol.* 32: 282–289.
6. Rauser *et al.* (2004) *Blood* 103: 3565–3572.

Antigen-specific T cells

IFN-γ-secreting antigen-specific T cells

Product	Regulatory status	Capacity (max.)	Availability	Order no.
CliniMACS Cytokine Capture System (IFN-gamma)	CE	1 × 10 ⁹ total cells	Europe, US ¹⁾	279-01
CliniMACS Cytokine Capture System (IFN-gamma) MicroBeads	For research use only, not for use in humans	1 × 10 ⁹ total cells	US, CAN, AUS	287-01

For availability in your country please contact your local representative.

1) In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

Description

Interferon-gamma (IFN-gamma) is a cytokine secreted by CD4⁺ and CD8⁺ memory and effector T cells upon antigenic restimulation.

The CliniMACS® Cytokine Capture System (IFN-gamma) is comprised of the CliniMACS IFN-gamma Catchmatrix Reagent, consisting of CD45 antibodies conjugated to IFN-gamma specific antibodies, and the CliniMACS IFN-gamma Enrichment Reagent. The CliniMACS IFN-gamma Enrichment Reagent consists of superparamagnetic iron dextran particles conjugated to IFN-gamma specific antibodies. The CliniMACS IFN-gamma Enrichment Reagent allows for magnetic selection of the IFN-gamma secreting cells.

One kit is suitable for processing up to 1 × 10⁹ total cells.

Please inquire about recommended device and accessories.

Applications

The CliniMACS® Cytokine Capture System (IFN-gamma) allows for the concomitant enrichment of viable antigen-specific CD4⁺ and CD8⁺ IFN-gamma secreting T cells.

Referenced literature

Antigen-specific CD4⁺ and CD8⁺ T cells can be enriched using the CliniMACS Cytokine Capture System (IFN-gamma)¹. T cells with specificity for CMV or Adenovirus could be expanded *in vitro*.²⁻³

The adoptive transfer of antigen-specific T cells may hold promise for the treatment of viral infections.⁴⁻⁵

References

1. Campbell (2003) *Methods* 31(2): 150–159.
2. Feuchtinger *et al.* (2004) *Exp. Hematol.* 32 (3): 282–289.
3. Rauser *et al.* (2004) *Blood* 103 (9): 3565–3572.
4. Feuchtinger *et al.* (2005) *Br. J. Haematol.* 128: 503–509.
5. Feuchtinger *et al.* (2006) *Br. J. Haematol.* 134: 64–76.

Related MACS® Products

CD4-FITC, human	130-080-501
CD4-APC, human	130-091-232
CD8-FITC, human	130-080-601
CD8-APC, human	130-091-076
Large-Scale IFN-γ Secretion Assay – Enrichment Kit, human	130-091-329
T cell Expansion Bag (6 ports)	130-091-459
T cell Expansion Bag (tube)	130-091-458
T cell Activation/Expansion Kit, human	130-091-441
Anti-IFN-γ-PE, human	130-091-653

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research only and not for therapeutic or diagnostic use.

Antigen-specific T cells

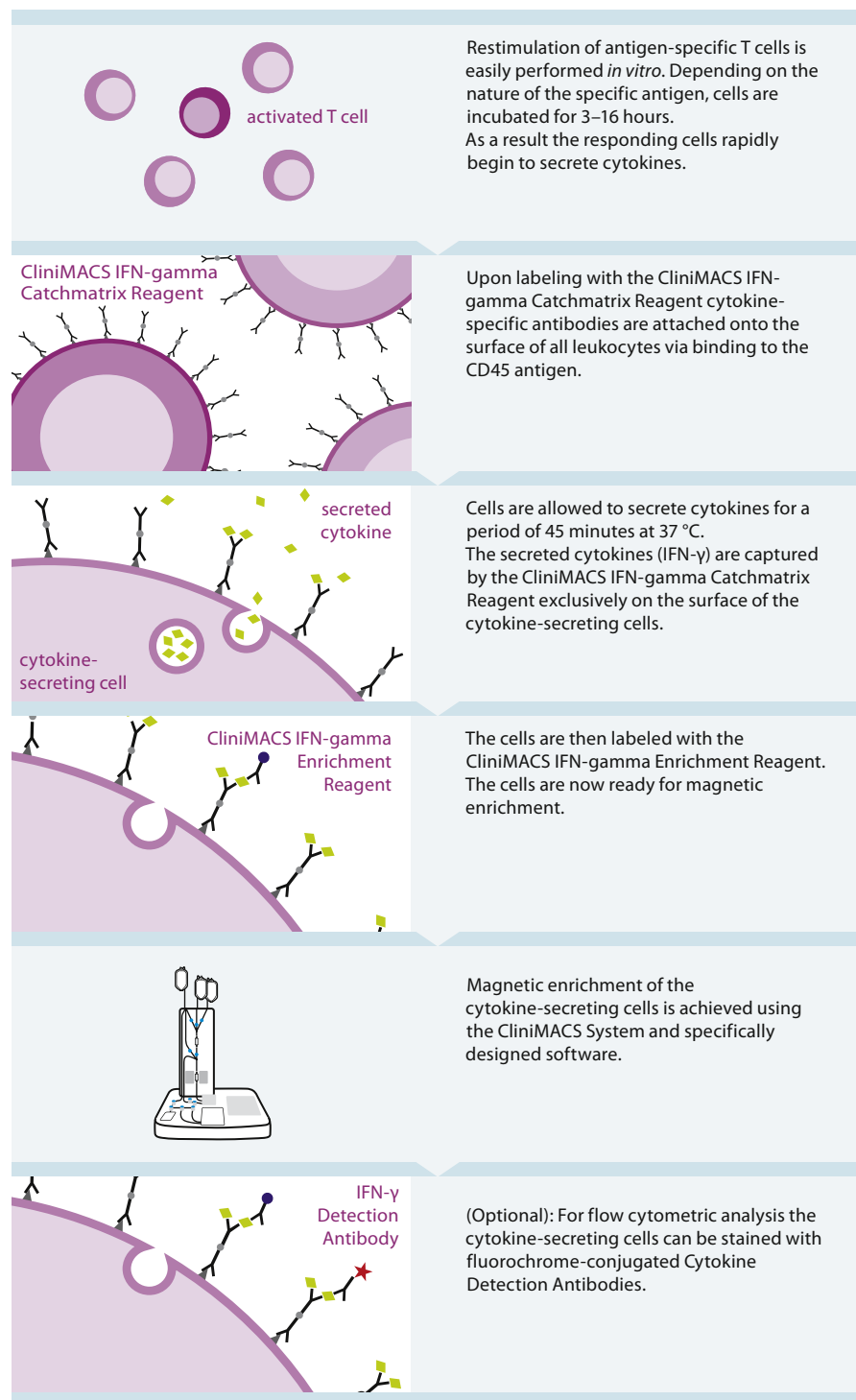


Figure 1: Principle of the CliniMACS® Cytokine Capture System (IFN-gamma).

NK cells

CliniMACS® Natural killer cell enrichment

Natural killer (NK) cells are known to be first in the line of defense against viral infections and solid tumors. On both malignant and virally infected cells a multiplicity of so called “danger signals” can be up-regulated. These signals can be specifically recognized by NK cells, enabling these cells to become activated and eventually lyse the potentially harmful target cell. It is the balance between stimulating and inhibiting signals which determines whether or not the NK cells will lyse their target cells¹. More recently, NK cells have been specifically associated with the induction of graft versus leukemia (GVL) effects in hematopoietic malignant diseases in murine models and some retrospective analyses in the human²⁻⁴. Based on these data, clinical studies⁵⁻⁷ were initiated to investigate the safety and feasibility of this approach in a haploidentical transplant setting for hematologic malignant disorders. The NK cells were enriched from leukapheresis products^{8,9} and given as donor lymphocyte infusions (DLI) at various time points post transplant. One important reason NK-DLI has been used instead of an unmanipulated DLI was the observation that NK cells may induce a GVL effect without inducing graft versus host disease (GVHD); the latter being a serious complication of unmanipulated DLIs in general. It has been theorized that a KIR-mismatch situation in the GVH direction may enable NK cells to lyse malignant target cells in order to exert GVL effects. This effect may be used to revert impending mixed chimerism or relapse in the above mentioned setting. Safety and feasibility of this approach was recently reviewed.¹⁰

References:

1. Kärre (2002) *Science* 295: 2029–2031.
2. Ruggeri *et al.* (2002) *Science* 295: 2097–2100.
3. Giebel *et al.* (2003) *Blood* 102: 814–819.
4. Elmaagacli *et al.* (2004) *Onkologie* 27, Suppl. 3: 0194.
5. Uharek *et al.* (2003) *Onkologie* 26, Suppl. 5: V374.
6. Passweg *et al.* (2003) *Onkologie* 26, Suppl. 5: P482.
7. Passweg *et al.* (2004) *Leukemia* 18 (11): 1835–1838.
8. Iyengar *et al.* (2003) *Cytotherapy* 5(6): 479–484.
9. Duwendag *et al.* (2004) 20th International NK cell Workshop, Noordwijkerhout, NL.
10. Passweg *et al.* (2005) *BMT* 35:637–643.

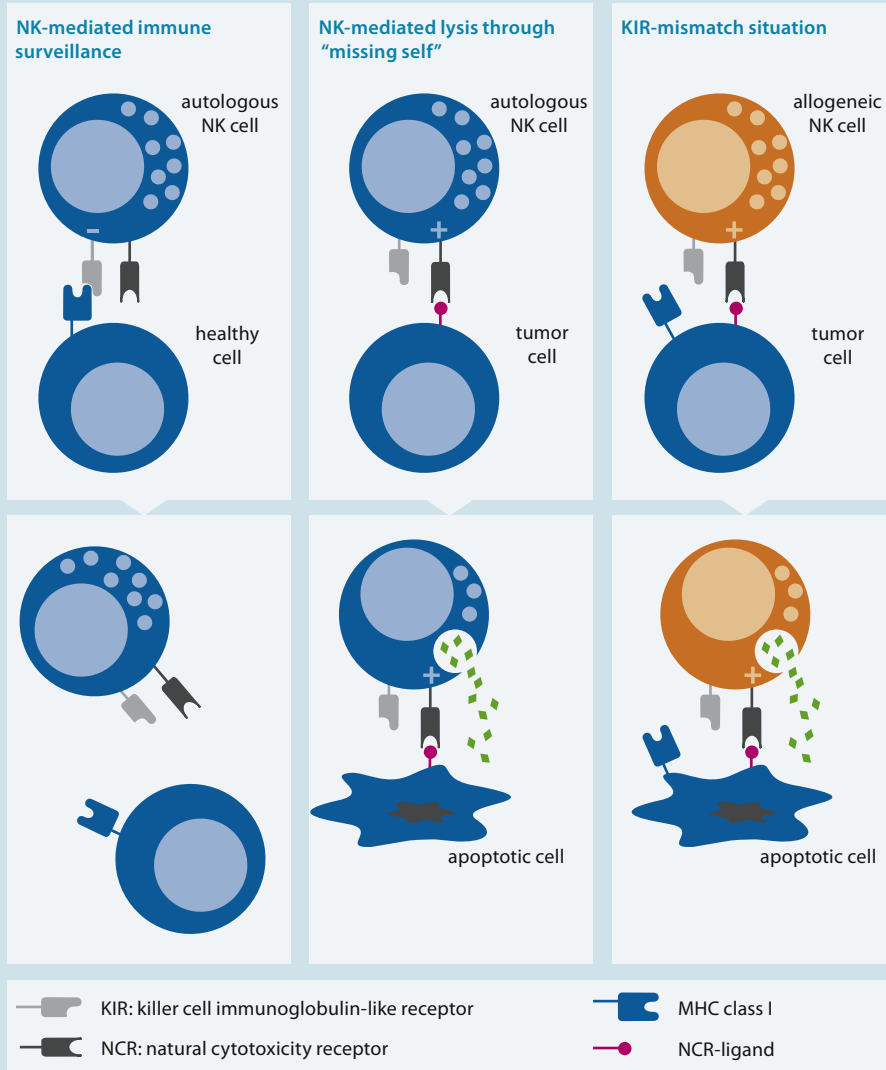


Figure 1: Defense mechanisms of natural killer cells.

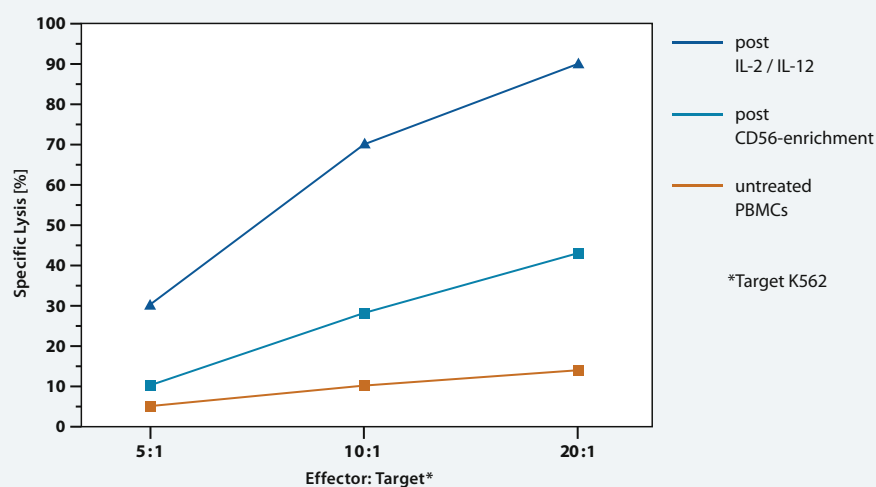


Figure 2: Activation of CD56⁺CD3⁻ NK cells upon treatment with IL-2 / IL-12. (Courtesy of Dr. Uharek, Berlin, Germany.)

CliniMACS® CD56 Cell Enrichment

Product	Regulatory status	Capacity (max.)	Availability	Order no.
CliniMACS CD56 Reagent	CE	10×10 ⁹ CD56 ⁺ cells from 40×10 ⁹ total cells	Europe, US ¹⁾	271-01
CliniMACS CD3/CD56 Reagent Combination	CE-marked components	15×10 ⁹ CD3 ⁺ cells and 10×10 ⁹ CD56 ⁺ cells from 40×10 ⁹ total cells	Europe	704-01
CliniMACS CD3/CD56 Complete Kit	CE-marked components	15×10 ⁹ CD3 ⁺ cells and 10×10 ⁹ CD56 ⁺ cells from 40×10 ⁹ total cells	Europe	740-04
CliniMACS CD56 MicroBeads	For research use only, not for use in humans	10×10 ⁹ CD56 ⁺ cells from 40×10 ⁹ total cells	US, CAN, AUS	194-01

For availability in your country please contact your local representative.

1) In the USA, the CliniMACS® System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

Description

The CD56 antigen is expressed on essentially all human NK cells, and its density on the cell membrane is increased after activation. The antigen is also present on a unique subset of CD3⁺ T cells (NKT cells) that mediate non-MHC-restricted cytotoxicity and on some neural tissues and tumors.

The CliniMACS® CD56 Products consist of superparamagnetic iron dextran particles directly conjugated to CD56 antibodies.

One vial contains 7.5 mL sterile, non-pyrogenic solution suitable for magnetic labeling of 10×10⁹ CD56⁺ cells from up to 40×10⁹ total cells.

Please inquire about recommended device and accessories.

Applications

CliniMACS® CD56 Products were developed for the enrichment of NK and NKT cells from human heterogeneous hematologic cell populations in combination with the CliniMACS® System.

In order to generate highly purified CD56⁺CD3⁺ NK cells, the CliniMACS CD56 Reagent is used in combination with the CliniMACS CD3 Reagent. In this approach, the cellular product is first depleted of CD3⁺ cells and then further enriched for CD56.^{1,2}

Referenced literature

NK cells have been used in haploidentical PBSC transplantation settings as donor lymphocyte infusions (DLIs) to explore the prevention of GVHD and early relapse of hematological malignant diseases (e.g. AML, CML).

Use of NK-DLIs for stabilizing an impending decreasing donor chimerism in a haploidentical transplantation setting has also been published.³⁻¹¹

References

1. Duwendag *et al.* (2004) 20th International NK Cell Workshop, Noordwijkerhout, NL.
2. Iyengar *et al.* (2003) *Cytotherapy*, 5(6): 479–484.
3. Uharek *et al.* (2002) *Onkologie*, 25, Suppl. 4: P844.
4. Uharek *et al.* (2003) *Onkologie*, 26, Suppl. 5: V374.
5. Passweg *et al.* (2003) *Onkologie*, 26, Suppl. 5: P482.
6. Leung *et al.* (2004) *J. Immunol.*, 172: 644–650.
7. Maillard *et al.* (2004) 8th International Symposium on DC, Brugge, Belgium.
8. Giebel *et al.* (2003) *Blood*, 102: 814–819.
9. Elmaagacli *et al.* (2004) *Onkologie* 27, Suppl. 3: O194.
10. Ruggeri *et al.* (2002) *Science*, 295: 2097–2100.
11. Farag *et al.* (2002) *Blood*, 100(6): 1935–1947.

NK cells

Kit components (not available in the US)

CliniMACS CD3/CD56 Reagent Combination (704-01):

- 1 vial CliniMACS CD3 Reagent
- 1 vial CliniMACS CD56 Reagent

CliniMACS CD3/CD56 Complete Kit (740-04):

- 1 vial CliniMACS CD3 Reagent
- 1 vial CliniMACS CD56 Reagent
- 1 CliniMACS Tubing Set
- 1 CliniMACS Depletion Tubing Set
- 2 units CliniMACS PBS/EDTA Buffer 3 × 1 L

Related CliniMACS® Products

CliniMACS CD3 MicroBeads	176-01
CliniMACS CD3 Reagent	273-01
CliniMACS CD3/CD19 Reagent Combination ¹⁾	704-00

1) Not available in the US.

Related MACS® Products

CD3-FITC, human	130-080-401
-----------------	-------------


The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research only and not for therapeutic or diagnostic use.

Single-step NK and NKT cell enrichment procedure

A) Before enrichment

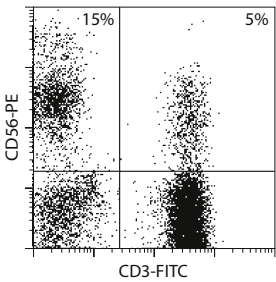

B) After enrichment


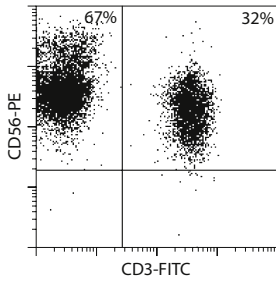



Figure 1: Flow cytometric analysis of CD56-enriched cells using the CliniMACS CD56 System.

Two-step NK cell enrichment procedure

A) Before enrichment

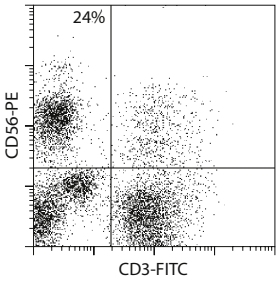

B) After CD3 depletion

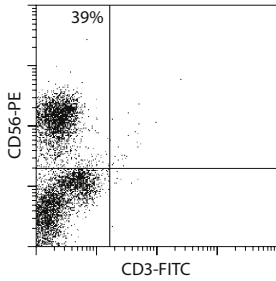

C) After CD3 depletion and CD56 enrichment

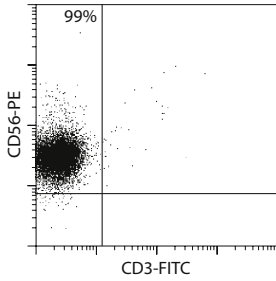


Figure 2: Flow cytometric analysis of CD3-depleted cells (B) followed by the analysis of CD56-enriched cells (C) using the CliniMACS CD3/CD56 System.

CliniMACS® CD19 Cell Depletion

Product	Regulatory status	Capacity (max.)	Availability	Order no.
CliniMACS CD19 Reagent	CE	5 × 10 ⁹ CD19 ⁺ cells from 40 × 10 ⁹ total cells	Europe, US ¹⁾	179-01
CliniMACS CD19 MicroBeads	For research use only, not for use in humans	5 × 10 ⁹ CD19 ⁺ cells from 40 × 10 ⁹ total cells	US, CAN, AUS	193-01

For availability in your country please contact your local representative.

1) In the USA, the CliniMACS® System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

Description

As a B cell lineage marker, CD19 is expressed from the early pro-B cell stage to the B cell lymphoblast stage; however, CD19 is down-regulated as B cells mature to plasma cells. The CD19 antigen is also expressed on most malignant B cells and on a subset of follicular dendritic cells.

The CliniMACS® CD19 Products consist of superparamagnetic iron dextran particles directly conjugated to CD19 antibodies.

One vial contains 7.5 mL sterile, non-pyrogenic solution suitable for magnetic labeling of up to 5 × 10⁹ CD19⁺ cells from up to 40 × 10⁹ total cells.

Please inquire about recommended device and accessories.

Applications


CliniMACS® CD19 Products were developed for depletion of CD19⁺ B cells from human heterogeneous hematologic cell populations in combination with the CliniMACS® System.

The depletion of CD19⁺ cells can be performed to remove unwanted malignant or potentially EBV-infected B cells from cellular products.¹

References

1. Barfield *et al.* (2004) Cytotherapy 6: 1–6.

A) Before depletion

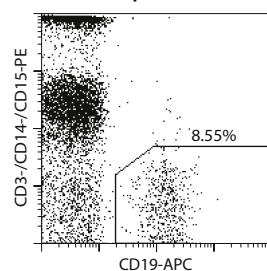

B) After depletion

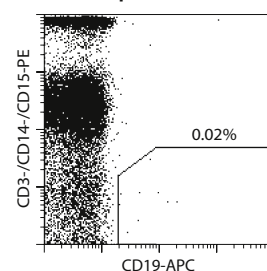


Figure 1: Flow cytometric analysis of CD19–depleted B cells from leukapheresis product using the CliniMACS CD19 System.

Related CliniMACS® Products

CliniMACS CD3/CD19 Reagent Combination ¹⁾	704-00
CliniMACS CD1c (BDCA-1)-Biotin	277-01

1) Not available in the US.

Related MACS® Products

CD20-FITC, human	130-091-108
CD3-PE, human	130-091-374
CD14-PE, human	130-091-242
CD15-PE, human	130-091-375

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research only and not for therapeutic or diagnostic use.

CliniMACS® Dendritic cell enrichment

Dendritic cells (DCs) have great potential as cellular vaccines for various diseases. Most clinical studies using DCs focus on malignant diseases such as solid tumors or hematologic malignancies. However, novel indications such as viral infections, solid organ transplantation, and autoimmune diseases are also being investigated.

DCs generated from monocytes after CD14 enrichment (MoDCs) have been used in many of these clinical trials and are well characterized.¹⁻⁵ *In vitro* experiments have demonstrated that MoDCs and DCs derived from CD34⁺ precursor cells are able to present antigens and efficiently induce a specific immune response dominated by effector T cells. However, in clinical studies using DCs, clinical responses have been lacking in most patients treated.

It is thought that other components of the immune system including NK cells, myeloid or plasmacytoid blood DCs (PDCs), as well as regulatory T cells may

be necessary for a clinical response. Recent findings support a role for regulatory T cells in limiting the tumor-specific immune response induced by DCs.⁶ It has been demonstrated that CD1c (BDCA-1)-positive myeloid DCs display a strong costimulatory and cross-presenting capacity.⁷

Plasmacytoid dendritic cells (PDCs) are the most potent Type I IFN-producing cells identified so far.⁸ Interferon-alpha-2a is used as pharmaceutical drug in indications such as malignant melanoma and other malignancies. Thus, the local application of PDCs could become a promising tool in enhanced vaccine strategies.

Recent developments will soon enable the use of blood myeloid DCs, which are CD1c (BDCA-1)-positive, or the use of CD304 (BDCA-4)-positive PDCs. The possibility to enrich both cell types will allow for new enhanced DC vaccine strategies. Because these DC subsets are derived directly from blood, no long culturing period – as for MoDC generation – is needed.

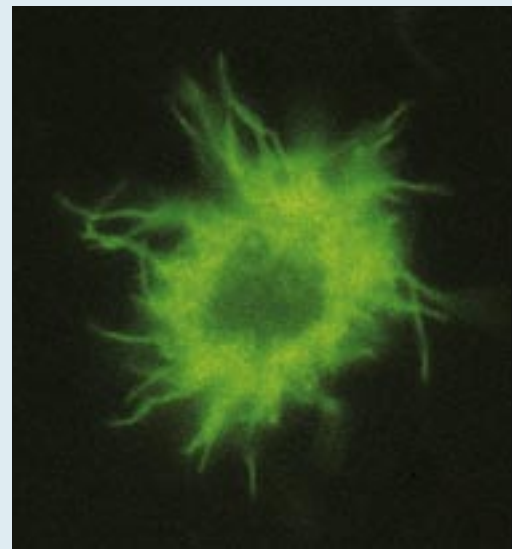


Figure 1: Mature blood dendritic cell stained with HLA-DR-FITC.

References:

1. Dietz *et al.* (2004) *Cytotherapy* 6(6): 563–570.
2. Curti *et al.* (2004) *Leukemia & Lymphoma* 45(7): 1419–1428.
3. Meyer-Wentrup *et al.* (2003) *J. Hematother. & Stem Cell Res.* 12: 289–299.
4. Sorg *et al.* (2003) *J. Immunol.* 26(4): 374–383.
5. Motta *et al.* (2003) *Br. J. Haematol.* 121: 240–250.
6. Prasad *et al.* (2005) *J. Immunol.* 174: 90–98.
7. Schnurr *et al.* (2005) *Blood* 105: 2465–2472.
8. Dzionek *et al.* (2002) *Hum. Immunol.* 63: 1133–1148.

CliniMACS® CD14 Cell Enrichment

Product	Regulatory status	Capacity (max.)	Availability	Order no.
CliniMACS CD14 Reagent	CE	4×10 ⁹ CD14 ⁺ cells from 20×10 ⁹ total cells	Europe, US ¹⁾	272-01
CliniMACS CD14 Complete Kit	CE-marked components	4×10 ⁹ CD14 ⁺ cells from 20×10 ⁹ total cells	Europe	281-01
CliniMACS CD14 MicroBeads	For research use only, not for use in humans	4×10 ⁹ CD14 ⁺ cells from 20×10 ⁹ total cells	US, CAN, AUS	191-01

For availability in your country please contact your local representative.

1) In the USA, the CliniMACS® System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

Description

The CD14 antigen belongs to the LPS receptor complex and monocytes strongly express the antigen.

The CliniMACS® CD14 Products consist of superparamagnetic iron dextran particles directly conjugated to CD14 antibodies.

One vial contains 7.5 mL sterile, non-pyrogenic solution suitable for magnetic labeling of up to 4×10⁹ CD14⁺ monocytes in 20×10⁹ total cells.


Please inquire about recommended device and accessories.

Applications

CliniMACS CD14 Products were developed for the enrichment of CD14⁺ cells from human heterogeneous hematologic cell populations in combination with the CliniMACS® System.

The enriched cells can be used for the subsequent generation of human monocyte-derived dendritic cells (MoDCs).¹

A) Before enrichment

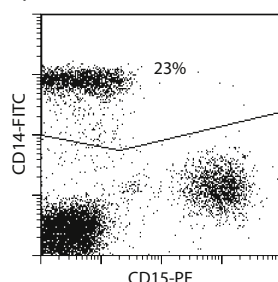

B) After enrichment

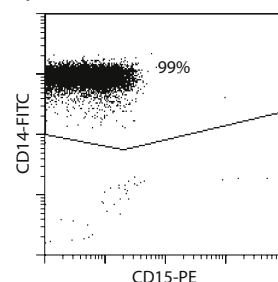


Figure 1: Flow cytometric analysis of CD14-enriched cells using the CliniMACS CD14 System.

Referenced literature

MoDCs are currently being used in clinical trials which mainly target malignant diseases such as solid tumors and hematologic malignancies.²⁻⁸

References

1. Campbell *et al.* (2005) *Methods in Molecular Medicine* 109 (5): 55–69.
2. Babatz *et al.* (2003) *J. Hematother. Stem Cell Res.* 12 (5): 515–523.
3. Motta *et al.* (2003) *Br. J. Hematol.* 121 (2): 240–250.
4. Meyer-Wentrup *et al.* (2003) *J. Hematother. Stem Cell Res.* 12 (3): 289–299.
5. Sorg *et al.* (2003) *J. Immunother.* 26 (4): 374–383.
6. Padley *et al.* (2001) *J. Hematother. Stem Cell Res.* 10: 427–429.
7. Dietz *et al.* (2000) *J. Hematother. Stem Cell Res.* 9 (1): 95–101.
8. Curti *et al.* (2004) *Leukemia & Lymphoma* 45 (7): 1419–1428.

Kit components (not available in the US)

CliniMACS CD14 Complete Kit (281-01):

- 1 vial CliniMACS CD14 Reagent
- 1 CliniMACS Tubing Set
- 3 × 1 L of CliniMACS PBS/EDTA Buffer

Related MACS® Products

CD14-FITC, human	130-080-701
CD15-FITC, human	130-081-101

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research only and not for therapeutic or diagnostic use.

CliniMACS® CD1c (BDCA-1) Cell Enrichment

Product	Regulatory status	Capacity (max.)	Availability	Order no.
CliniMACS CD1c (BDCA-1)-Biotin	CE	0.2×10^9 CD1c (BDCA-1) ⁺ cells from 40×10^9 total cells	Europe, US ¹⁾	NEW 277-01
CliniMACS CD1c (BDCA-1)/CD19 Complete Kit	CE-marked components	0.2×10^9 CD1c (BDCA-1) ⁺ cells from 40×10^9 total cells	Europe	NEW 740-03
CliniMACS CD1c (BDCA-1)-Biotin for Research Use	For research use only, not for use in humans	0.2×10^9 CD1c (BDCA-1) ⁺ cells from 40×10^9 total cells	US, CAN, AUS	255-01

For availability in your country please contact your local representative.

1) In the USA, the CliniMACS® System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

Description

The CD1c (BDCA-1) antigen is expressed on dendritic cells which are CD11c^{high}CD123^{low} and represent a major subset of myeloid dendritic cells in human peripheral blood. Also, the antigen is expressed on a subpopulation of CD19⁺ B cells.

The CliniMACS® CD1c (BDCA-1)-Biotin Products consist of biotin directly conjugated to CD1c (BDCA-1) antibodies.

One vial contains 5 mL sterile, non-pyrogenic solution suitable for labeling of up to 0.2×10^9 CD1c (BDCA-1)⁺ cells from up to 40×10^9 total cells.

Please inquire about recommended device and accessories.

Applications

CliniMACS CD1c (BDCA-1)-Biotin Products were developed for the enrichment of CD1c (BDCA-1)⁺ myeloid dendritic cells from human heterogeneous hematologic cell populations in combination with the CliniMACS® System.

After labeling of target cells with the CliniMACS CD1c (BDCA-1)-Biotin, the CliniMACS Anti-Biotin Reagent is used for the magnetic enrichment of CD1c (BDCA-1)⁺ cells from the cellular product.

A CD19⁺ B cell depletion is recommended prior to the enrichment of CD1c (BDCA-1)⁺ cells.

Flow cytometric analysis of CD1c (BDCA-1)⁺ cell enrichment

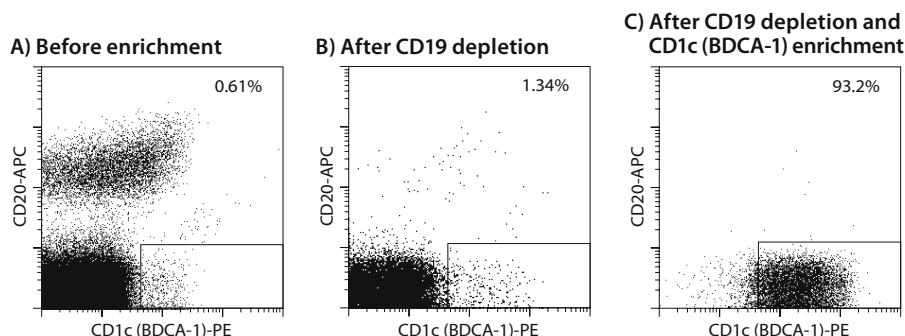


Figure 1: Flow cytometric analysis of CD19-depleted cells (B) followed by the analysis of CD1c (BDCA-1)-enriched cells (C) using the CliniMACS CD19- and CliniMACS CD1c (BDCA-1)-Biotin System in combination.

Kit components (not available in the US)

CliniMACS CD1c (BDCA-1)/CD19 Complete Kit (740-03):

- 1 vial CliniMACS CD1c (Anti-BDCA-1)-Biotin
- 1 vial CliniMACS CD19 Reagent
- 1 vial CliniMACS Anti-Biotin Reagent
- 1 CliniMACS Tubing Set
- 1 CliniMACS Tubing Set LS
- 3 units 3x1L CliniMACS PBS/EDTA Buffer

References

1. Dzionek A. *et al.* (2000) J. Immunol. 165: 6037–6046.
2. Dzionek A. *et al.* (2002) Hum Immunol. 63 (12): 1133–1148.
3. Piechaczek C *et al.* (2002) 7th International Symposium on Dendritic Cells, Bamberg, Germany: P197.
4. Schnurr M. *et al.* (2005) Blood 105(6): 2465–2472.
5. Davis I.D. *et al.* (2005) Proc. Amer. Assoc. Cancer Res. 46: Abstr. no. 3466.

Related CliniMACS® Products

CliniMACS Anti-Biotin MicroBeads	192-01
CliniMACS Anti-Biotin Reagent	173-01
CliniMACS CD19 MicroBeads	193-01
CliniMACS CD19 Reagent	179-01

Related MACS® Products

Blood Dendritic Cell Enumeration Kit, human	130-091-324
CD1c (BDCA-1)-PE, human	130-090-508

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research only and not for therapeutic or diagnostic use.

CliniMACS® CD304 (BDCA-4) Cell Enrichment

Product	Regulatory status	Capacity (max.)	Availability	Order no.
CliniMACS CD304 (BDCA-4) MicroBeads	For research use only, not for use in humans	0.2 × 10 ⁹ CD304 (BDCA-4) ⁺ cells from 40 × 10 ⁹ total cells	Worldwide	291-01

For availability in your country please contact your local representative.

Description

In peripheral blood, CD304 (BDCA-4, Neuropilin-1) is expressed on human plasmacytoid (lymphoid) dendritic cells, which are characterized as being CD303 (BDCA-2)⁺, CD4⁺, Lin⁻, CD11c⁻, CD123^{bright}, CD45RA⁺, CD2⁻. These cells do not express myeloid lineage markers such as CD13 or CD33^{1,2}.

The CliniMACS® CD304 (BDCA-4) MicroBeads consist of superparamagnetic iron dextran particles directly conjugated to CD304 (BDCA-4) antibodies.

One vial contains 7.5 mL sterile, non-pyrogenic solution suitable for magnetic labeling of up to 0.2 × 10⁹ CD304 (BDCA-4)⁺ cells out of 40 × 10⁹ total cells.

Please inquire about recommended device and accessories.


Applications

CliniMACS® CD304 (BDCA-4) MicroBeads were developed for enrichment of CD304 (BDCA-4)⁺ plasmacytoid dendritic cells from human heterogeneous hematologic cell populations in combination with the CliniMACS® System.

References

1. Dzionek *et al.* (2000) *J. of Immunology* 165: 6037–6046. [898]
2. Dzionek *et al.* (2002) *Hum Immunol.* 63 (12): 1133–1148. [2423]

A) Before enrichment

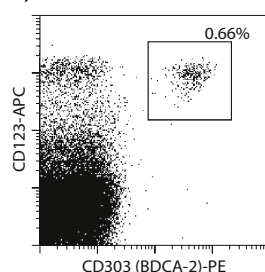

B) After enrichment

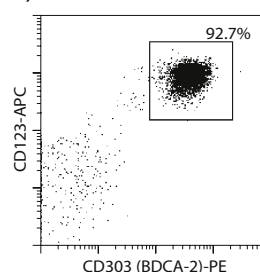


Figure 1: Flow cytometric analysis of CD304 (BDCA-4)-enriched cells from leukapheresis harvest using the CliniMACS CD304 (BDCA-4) System.

Related MACS® Products

CD303 (BDCA-2)-FITC, human	130-090-510
CD303 (BDCA-2)-PE, human	130-090-511
CD123-APC, human	130-090-901
CD45-FITC, human	130-080-202
CD45-PE, human	130-080-201
Blood Dendritic Cell Enumeration Kit, human	130-091-324

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research only and not for therapeutic or diagnostic use.

CliniMACS® Anti-Biotin

Product	Regulatory status	Capacity	Availability	Order no.
CliniMACS Anti-Biotin Reagent	CE	40×10 ⁹ total cells	Europe, US ¹⁾	173-01
CliniMACS Anti-Biotin MicroBeads	For research use only, not for use in humans	40×10 ⁹ total cells	US, CAN, AUS	192-01

For availability in your country please contact your local representative.

1) In the USA, the CliniMACS® System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

Description

The CliniMACS® Anti-Biotin Products consist of superparamagnetic iron dextran particles directly conjugated to Biotin-specific antibodies.

One vial contains 7.5 mL sterile, non-pyrogenic solution suitable for magnetic labeling of target cells from up to 40×10⁹ total cells. Please inquire about recommended device and accessories.

Applications

The CliniMACS® Anti-Biotin Products were developed for the enrichment or depletion of target cells labeled with a biotinylated primary antibody or ligand.

Related CliniMACS® Products

CliniMACS CD1c (BDCA-1)-Biotin 277-01

Related MACS® Products

Anti-Biotin-FITC	130-090-857
Anti-Biotin-PE	130-090-756
Anti-Biotin-APC	130-090-856

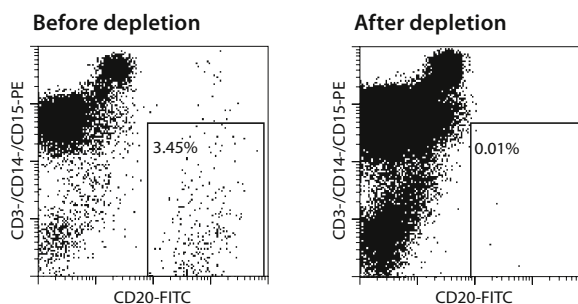


Figure 1: Flow cytometric analysis of CD19-depleted cells using the CliniMACS Anti-Biotin System in combination with CD19-biotinylated antibody.

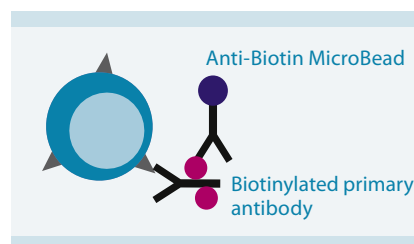


Figure 2: Labeling principle using CliniMACS Anti-Biotin Products.

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research only and not for therapeutic or diagnostic use.

CliniMACS® Biotinylation Kit 500

Product	Regulatory status	Capacity	Availability	Order no.
CliniMACS Biotinylation Kit 500 ¹⁾	CE	Biotinylation of 500 µg monoclonal antibody	Europe, US ²⁾ NEW	701-06
CliniMACS Biotinylation Kit 500 for Research Use	For research use only, not for use in humans	Biotinylation of 500 µg monoclonal antibody	US, CAN, AUS	257-01

For availability in your country please contact your local representative.

1) Outside Europe the CliniMACS® Biotinylation Kit 500 is for research use only.

2) In the USA, the CliniMACS® System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

Description/Applications

The CliniMACS® Biotinylation Kit 500 Products have been developed for the biotinylation of monoclonal antibodies for further magnetic cell enrichment from human heterogeneous hematologic cell populations using the CliniMACS System.

One kit is sufficient for biotinylation 500 µg monoclonal antibody.

Please check in advance whether the antibody is suitable for biotinylation, magnetic labeling, and cell separation using the CliniMACS® System.

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).

CliniMACS® MicroBeads are for research use only and not for use in humans.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research only and not for therapeutic or diagnostic use.

CliniMACS® PBS/EDTA Buffer				
Product	Content	Regulatory status	Availability	Order no.
CliniMACS PBS/EDTA Buffer	3 × 1L plastic bags	CE	Europe, CAN, AUS, US ¹⁾	700-25
For availability in your country please contact your local representative.				
1) In the USA, the CliniMACS® System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE).				

Description

The CliniMACS® PBS/EDTA Buffer consists of phosphate buffered saline, pH 7.2, supplemented with 1 mM EDTA. The buffer is supplied in 1 L plastic bags as a sterile, non-pyrogenic solution.

The CliniMACS® System components: Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS System components are available as CE-marked medical devices. In the USA, the CliniMACS System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE). CliniMACS® MicroBeads are for research use only and not for use in humans.



Figure 1: CliniMACS PBS/EDTA Buffer.

Application

CliniMACS PBS/EDTA Buffer was developed for cell preparation and separation of magnetically labeled cells with the CliniMACS System. CliniMACS PBS/EDTA Buffer is not intended for direct infusion into humans.

CliniMACS® PBS/EDTA Buffer for Research Use				
Product	Content	Regulatory status	Availability	Order no.
CliniMACS PBS/EDTA Buffer for Research Use	3 × 1L plastic bags	For research use only, not for use in humans ¹⁾	US	705-25
For availability in your country please contact your local representative.				
1) Manufactured under an ISO 9001 certified quality system.				

Description

The CliniMACS® PBS/EDTA Buffer consists of phosphate buffered saline, pH 7.2, supplemented with 1 mM EDTA. The buffer is supplied in 1 L plastic bags as a sterile, non-pyrogenic solution.

Application

CliniMACS PBS/EDTA Buffer was developed for cell preparation and separation of magnetically labeled cells with the CliniMACS® System.



Figure 1: CliniMACS PBS/EDTA Buffer for research use.

Cell culture products

The technical features of Miltenyi Biotec cell culture bags make them valuable tools for investigators interesting in culturing cells.

The bags are gas permeable, transparent for microscopy and allow for a closed culture system resulting in reduced risk of contamination.

The bags are individually packed, sterile and tested for endotoxins.

They come in single-tube format or with six needleless access connectors fitting standard laboratory requirement.

The single-tube can be used for sterile docking (according to the instructions for use of the tube welding device).


Transfer devices with a male luer lock connector (e.g. tubing or syringe) can be attached directly to the needleless access connectors.

Cell Differentiation Bags are available for three different culture volumes: 100 mL, 250 mL and 500 mL.

Cell Expansion Bags are compartmentalized with easy to open seals for expandable culture volumes as cell numbers increase during culture period. There is no need for changing culture vessel.

The three compartments allow for expandable volumes from 8 mL up to 100 mL.

1.

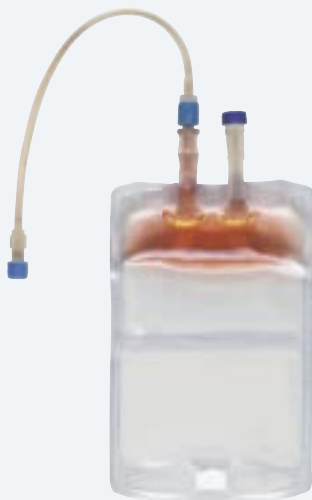

2.



3.



Cell Expansion Bag (tube) with incremental culture volumes: As culture volume increases, seals can easily be opened to give larger culture volumes. No transfer to other bags is necessary.

Cell Expansion Bags						
Product	Content	Regulatory status	Capacity	Availability	Order no.	
Cell Expansion Bag (tube)	5 Bags	CE ¹⁾	up to 100 mL	Europe	200-074-301	
Cell Expansion Bag (6 ports)	5 Bags	CE ¹⁾	up to 100 mL	Europe	200-074-302	
Cell Expansion Bag (tube)	5 Bags	For research use only ²⁾	up to 100 mL	US, CAN, AUS	NEW	130-074-351
Cell Expansion Bag (6 ports)	5 Bags	For research use only ²⁾	up to 100 mL	US, CAN, AUS	NEW	130-074-352

For availability in your country please contact your local representative.

1) The Cell Expansion Bags are manufactured by Miltenyi Biotec GmbH and controlled under an ISO 13485 certified quality system. These products are available in Europe as CE-marked medical devices.

2) Manufactured by Miltenyi Biotec GmbH under an ISO 9001 certified quality system.

Description

Cell Expansion Bags are gas-permeable and transparent for microscopy.

They have three compartments to allow for expandable volumes of up to 100 mL.
The bag are individually packed, sterile and tested for endotoxins.

Applications

Cell Expansion Bags are intended for the *in vitro* cultivation and expansion of human cells from heterogeneous haematologic cell populations (e.g. for the expansion of antigen-specific T cells).

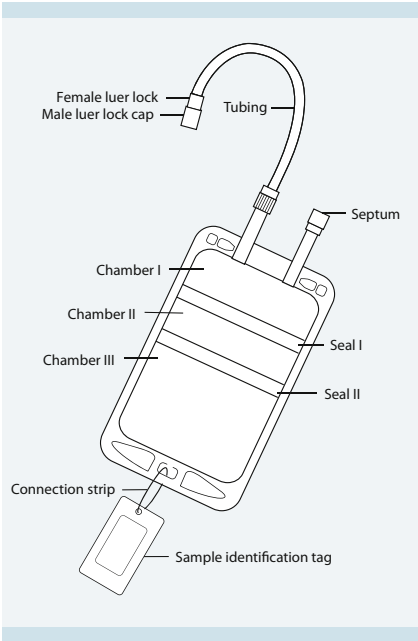


Figure 1: Cell Expansion Bag (tube) with two seals for expandable culture volumes.

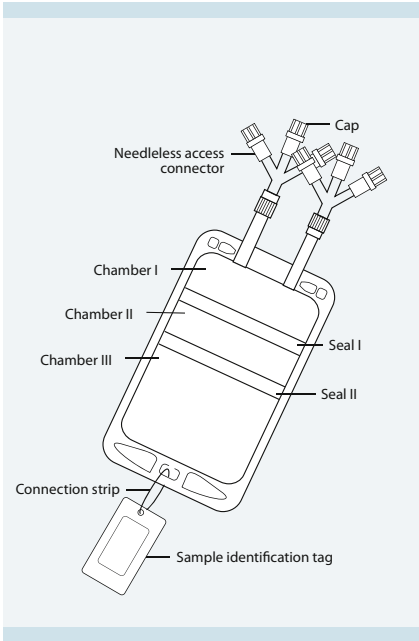


Figure 2: Cell Expansion Bag (6 ports) with two seals for expandable culture volumes and six needleless access connectors.

Accessories

Cell Differentiation Bags

Product	Content	Regulatory status	Capacity	Availability	Order no.
Cell Differentiation Bag (tube) – 100	5 Bags	CE ¹⁾	100 mL	Europe	200-074-101
Cell Differentiation Bag (tube) – 250	5 Bags	CE ¹⁾	250 mL	Europe	200-074-102
Cell Differentiation Bag (tube) – 500	2 Bags	CE ¹⁾	500 mL	Europe	200-074-103
Cell Differentiation Bag (6 ports) – 100	5 Bags	CE ¹⁾	100 mL	Europe	200-074-201
Cell Differentiation Bag (6 ports) – 250	5 Bags	CE ¹⁾	250 mL	Europe	200-074-202
Cell Differentiation Bag (6 ports) – 500	2 Bags	CE ¹⁾	500 mL	Europe	200-074-203
Cell Differentiation Bag (tube) – 100	5 Bags	For research use only ²⁾	100 mL	US, CAN, AUS NEW	130-074-151
Cell Differentiation Bag (tube) – 250	5 Bags	For research use only ²⁾	250 mL	US, CAN, AUS NEW	130-074-152
Cell Differentiation Bag (tube) – 500	2 Bags	For research use only ²⁾	500 mL	US, CAN, AUS NEW	130-074-153
Cell Differentiation Bag (6 ports) – 100	5 Bags	For research use only ²⁾	100 mL	US, CAN, AUS NEW	130-074-251
Cell Differentiation Bag (6 ports) – 250	5 Bags	For research use only ²⁾	250 mL	US, CAN, AUS NEW	130-074-252
Cell Differentiation Bag (6 ports) – 500	2 Bags	For research use only ²⁾	500 mL	US, CAN, AUS NEW	130-074-253

For availability in your country please contact your local representative.

1) The Cell Differentiation Bags are manufactured and controlled under an ISO 13485 certified quality system. These products are available in Europe as CE-marked medical devices.

2) Manufactured by Miltenyi Biotec GmbH under an ISO 9001 certified quality system.

Description

Cell Differentiation Bags are gas-permeable and transparent for microscopy.

They are available for three different culture volumes: 100 mL, 250 mL and 500 mL.

The bags are individually packed, sterile and tested for endotoxins.

Applications

Cell Differentiation Bags are intended for the *in vitro* cultivation and expansion of cells from heterogeneous haematologic cell populations (e.g. for the generation of dendritic cells from CD14-enriched monocytes).

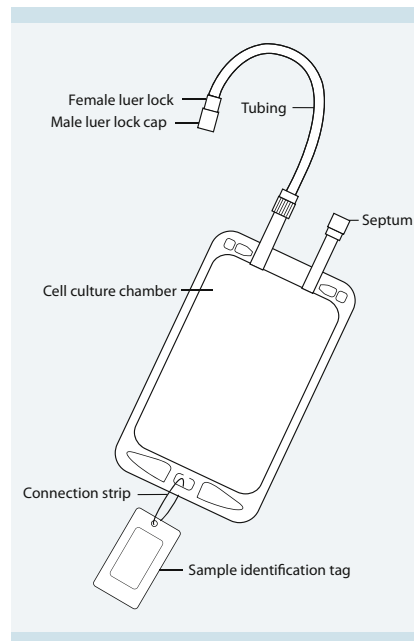


Figure 1: Cell Differentiation Bag (tube).

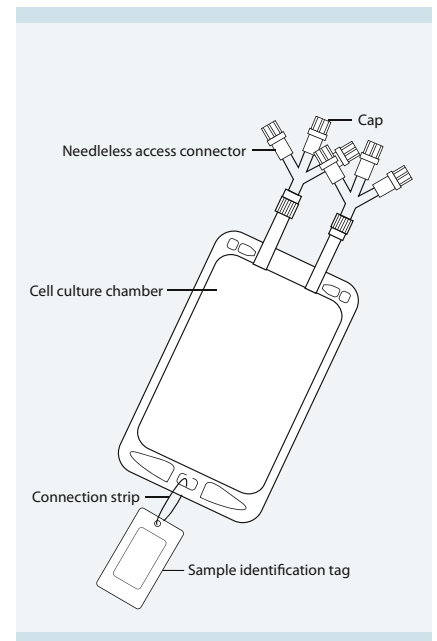


Figure 2: Cell Differentiation Bag (6 ports) with six needleless access connectors.

Pre-System Filter

Product	Content	Regulatory status	Availability	Order no.
Pre-System Filter	5 Filters	CE	Europe	181-01

For availability in your country please contact your local representative.

Description

Pre-System Filter for cell separation (Blood transfusion filter, 40 micron pore width). A Pre-System Filter should be used with all CliniMACS® Tubing Sets in order to retain cell aggregates.

Sterilized, non-pyrogenic fluid pathway, single-packed.

Please inquire about recommended device and accessories.



Figure 1: The Pre-System Filter.

Luer/Spike Interconnector

Product	Content	Regulatory status	Availability	Order no.
Luer/Spike Interconnector	5 Interconnectors	CE	Europe	187-01

For availability in your country please contact your local representative.

Description

Adapter with spike and female luer connector, suitable for the attachment of transfer bags to these CliniMACS® Tubing Sets.

Sterilized, non-pyrogenic, single-packed.

Please inquire about recommended device and accessories.



Figure 1: The Luer/Spike Interconnector.

Sampling Site Coupler

Product	Content	Regulatory status	Availability	Order no.
Sampling Site Coupler	5 Couplers	CE	Europe	189-01

For availability in your country please contact your local representative.

Description

Injection port, suitable for cell labeling procedures (e.g. for taking samples). Sampling Site Coupler is suitable for centrifugation (connected to transfer bags).

Sterilized, single-packed.

Please inquire about recommended device and accessories.



Figure 1: The Sampling Site Coupler.

Transfer Set Coupler/Needle

Product	Content	Regulatory status	Availability	Order no.
Transfer Set Coupler/Needle	5 Transfer Sets	CE	Europe	185-01

For availability in your country please contact your local representative.

Description

Plasma transfer set with coupler and puncture needle.

This transfer set can be used for connecting the CliniMACS® PBS/EDTA Buffer bag to the cell preparation bag.

Sterilized, non-pyrogenic, single-packed.

Please inquire about recommended device and accessories.



Figure 1: The Transfer Set Coupler/Needle.

Transfer Set Coupler/Coupler

Product	Content	Regulatory status	Availability	Order no.
Transfer Set Coupler/Coupler	5 Transfer Sets	CE	Europe	186-01

For availability in your country please contact your local representative.

Description

Plasma transfer set with two couplers.

This transfer set can be used for connecting the CliniMACS® PBS/EDTA Buffer bag to the cell preparation bag or to connect two buffer bags.

Sterilized, non-pyrogenic, single-packed.

Please inquire about recommended device and accessories.



Figure 1: The Transfer Set Coupler/Coupler.

Transfer Set Needle/Needle

Product	Content	Regulatory status	Availability	Order no.
Transfer Set Needle/Needle	5 Transfer Sets	CE	Europe	188-01

For availability in your country please contact your local representative.

Description

Plasma transfer set with two needles.

This transfer set can be used for connecting the CliniMACS® PBS/EDTA Buffer Bag to the Cell Preparation Bag.

Sterilized, non-pyrogenic, single-packed.

Please inquire about recommended device and accessories.



Figure 1: The Transfer Set Needle/Needle.

Transfer Bags 150 mL, 600 mL, and 1000 mL

Product	Content	Regulatory status	Availability	Order no.
Transfer Bag 150 mL	5 Bags	CE	Europe	183-01
Transfer Bag 600 mL	5 Bags	CE	Europe	190-01
Transfer Bag 1000 mL	5 Bags	CE	Europe	180-01

For availability in your country please contact your local representative.

Description

Transfer Bag 150 mL

Suitable for collecting the positive fraction during a CliniMACS® Separation (in combination with a Luer/Spike Interconnector).

Sterilized, non-pyrogenic, single-packed.

Transfer Bag 600 mL

Suitable for cell labeling procedures.

Sterilized, non-pyrogenic, single-packed.

Transfer Bag 1000 mL

Suitable for collecting fractions during a CliniMACS® Separation (in combination with a Luer/Spike Interconnector).

Sterilized, non-pyrogenic, single-packed.

Please inquire about recommended device and accessories.



Figure 1: The Transfer Bag 150 mL.

Transfer Bag 600 mL with 8 couplers

Product	Content	Regulatory status	Availability	Order no.
Transfer Bag 600 mL with 8 couplers	5 Bags	CE	Europe	184-01

For availability in your country please contact your local representative.

Description

The Transfer Bag with eight couplers ("Octopus Bag") is suitable for cell labeling procedures.

This Transfer Bag enables the operator to perform the sample preparation procedure in a closed system.

Sterilized, non-pyrogenic, single-packed.

Please inquire about recommended device and accessories.



Figure 1: The Transfer Bag 600 mL with 8 couplers ("Octopus bag").

Distributors: products for clinical use

Argentina	Lab Systems S.A.	Quesada 5916 CP 1431 Buenos Aires Ciudad Autónoma de Buenos Aires , Argentina	Phone +54 11 4572 8458 Fax +54 11 4572 8458	direccion@lab-systems.com.ar www.lab-systems.com.ar
Australia	Miltenyi Biotec Australia Pty. Ltd.	Unit 16 A 2 Eden Park Drive North Ryde, NSW 2113, Australia	Phone +61 02 8877 7400 Fax +61 02 9889 5044	macs@miltenyibiotec.com.au www.miltenyibiotec.com
Austria	Miltenyi Biotec GmbH	Friedrich-Ebert-Straße 68 51429 Bergisch Gladbach Germany	Phone +49 2204 8306-0 Fax +49 2204 85197	macs@miltenyibiotec.de www.miltenyibiotec.com
Bahrain	Al Yamama	Al Saif Medical Centre Ibn Khaldoun Street Amman, Jordan	Phone +962 6 4655 166 Fax +962 6 4655 167	modhafar@alyamamagroup.com www.alyamamagroup.com
Belarus	Medbioscience	Nezavisimosti av., 107-11 220114 Minsk, Belarus	Phone +375 172 67 00 25 Fax +375 172 67 74 32	bioscience@open.by
Belgium	Miltenyi Biotec B.V.	Postbus 85183 3508 AD Utrecht The Netherlands	Phone (Belgium) 0800 94016 Fax (Belgium) 0800 99626	macs@miltenyibiotec.nl www.miltenyibiotec.com
Brazil	Ambriex S/A	Rua-Tupi, 535 - Santa Cecília 01233-907 Sao Paulo, Brazil	Phone +55 11 3665 1090 Fax +55 11 3665 1096	bio1@ambriex.com.br www.ambriex.com.br
Bulgaria	Epsilon & Epsilon Medical S.A.	Apollon Tower, 15 A3 64 L. Riankour str. 11523 Athens, Greece	Phone +30 210 6996 191 Fax +30 210 6920 210	info@eemed.gr www.eemed.gr
Canada	Miltenyi Biotec Inc.	12740 Earhart Avenue Auburn, CA 95602, USA	Phone +1 530 888 88 71 Fax +1 530 888 89 25	macs@miltenyibiotec.com www.miltenyibiotec.com
Chile	UB Ursula Biggemann y Cía Ltda.	Vitacura 2909 Of. 1007 Santiago – Las Condes, Chile	Phone +56 2 334 9282 Fax +56 2 334 1136	ursulab@entelchile.net www.ub.co.cl
China	Miltenyi Biotec Trading (Shanghai) Co., Ltd., Shanghai Office	Rm. 2301, No. 319 Xianxia Rd. Fareast International Plaza A Shanghai 200051, P. R. China	Phone +86 21 6235 1005 Fax +86 21 6235 0953	macs@miltenyibiotec.com.cn www.miltenyibiotec.com
	Miltenyi Biotec Trading (Shanghai) Co., Ltd., Beijing Office	Unit 916, Silver Tower, #2 Dong San Huan North Road Chao Yang District Beijing 100027, P. R. China	Phone +86 10 6410 7101 Fax +86 10 6410 7102	macs@miltenyibiotec.com.cn www.miltenyibiotec.com
Colombia	Equimed Ltda.	Calle 37 #16-51 , Bogotá DC Colombia	Phone +57 1 285 5053 Fax +57 1 285 1538	equimedcomercial@etb.net.co
Croatia	MEDIVA d.o.o.	Savska cesta 141 10000 Zagreb, Croatia	Phone +385 1 6191 580 Fax +385 1 6191 608	damir.jambrosic@mediva.hr www.mediva.hr
Cyprus	Epsilon & Epsilon Medical S.A.	Apollon Tower, 15 A3 64 L. Riankour str. 11523 Athens, Greece	Phone +30 210 6996 191 Fax +30 210 6920 210	info@eemed.gr www.eemed.gr
Czech Republic	BIOHEM spol.s.r.o.	Zlátovska 2211 911 01 Trenčín, Slovakia	Phone +421 32 6524 998 Fax +421 32 6525 792	biohem@biohem.sk www.biohem.sk
Ecuador	Frisonex	C. La Tierra #388/E8-11B y Av. de los Shyris 17-17-1413 Casilla, Ecuador	Phone +593 22 267 554 Fax +593 22 433 550	frisonex@frisonex.com
Egypt	Al Yamama	Al Saif Medical Centre Ibn Khaldoun Street Amman, Jordan	Phone +962 6 4655 166 Fax +962 6 4655 167	modhafar@alyamamagroup.com www.alyamamagroup.com

Distributors: products for clinical use

Estonia	AS Oriola	Department of Laboratory Supplies & Equipment Kungla Street 2 76505 Saue, Estonia	Phone +372 65 15 159 Fax +372 65 15 111	kaari.saarma@oriola.com www.oriola.ee
France	Miltenyi Biotec SAS	10 rue Mercœur 75011 Paris, France	Phone +33 1 56 98 16 16 Fax +33 1 56 98 16 17	macs@miltenyibiotec.fr www.miltenyibiotec.com
Germany	Miltenyi Biotec GmbH	Friedrich-Ebert-Straße 68 51429 Bergisch Gladbach Germany	Phone +49 2204 8306-0 Fax +49 2204 85197	macs@miltenyibiotec.de www.miltenyibiotec.com
Greece	Epsilon & Epsilon Medical S.A.	Apollon Tower, 15 A3 64 L. Riankour str. 11523 Athens, Greece	Phone +30 210 6996 191 Fax +30 210 6920 210	info@eemed.gr www.eemed.gr
Hong Kong	Miltenyi Biotec Trading (Shanghai) Co., Ltd., Shanghai Office	Rm. 2301, No. 319 Xianxia Rd. Fareast International Plaza A Shanghai 200051, P. R. China	Phone +86 21 6235 1005 Fax +86 21 6235 0953	macs@miltenyibiotec.com.cn www.miltenyibiotec.com
Hungary	Frank Diagnosztika Kft.	Dereglye u. 2 1036 Budapest, Hungary	Phone +361 388 3114 Fax +361 368 5721	frankdiagn@frank-diagn.hu www.frank-diagn.hu
India	Medicons	GH-8/505 Paschim Vihar New Delhi-87, India	Phone +91 11 6543 8094 Fax +91 11 2528 7644	medicons@mediconsinc.com www.mediconsinc.com
Indonesia	PT. TRANSMEDIC INDONESIA	Graha Atrium Lt. 6 Suite 6.02A Jalan Senen Raya No. 135 Jakarta Pusat 10410, Indonesia	Phone +62 21 3509 544 Fax +62 21 3509 545	www.transmedic.com.sg doni@transmedic-ind.com
Israel	Y.A. Almog Diagnostic & Medical Equipment Ltd.	North industrial area 4 60850 Shoham, Israel	Phone +972 3 977 33 90 Fax +972 3 977 33 91	info@almog.co.il www.almog.co.il
Italy	Miltenyi Biotec S.r.l.	via Persicetana, 2/D 40012 Calderara di Reno (Bologna), Italy	Phone +39 051 64 60 411 Fax +39 051 64 60 499	macs@miltenyibiotec.it www.miltenyibiotec.com
Japan	Miltenyi Biotec K.K.	Nittsu-Eitai Building 5F 16-10 Fuyuki, Koto-ku 135-0041 Tokyo, Japan	Phone +81 3 5646 8910 Fax +81 3 5646 8911	macs@miltenyibiotec.jp www.miltenyibiotec.co.jp
Jordan	Al Yamama	Al Saif Medical Centre Ibn Khaldoun Street Amman, Jordan	Phone +962 6 4655 166 Fax +962 6 4655 167	modhafar@alyamamagroup.com www.alyamamagroup.com
Kuwait	TRI ALPHA CO.	Al-Awadi Towers, Tower No. 3, 6th. Floor, Ahmend Al-Jaber Street, 72263 Da. A. Al-Salem, Kuwait	Phone +965 24 3 1344 Fax +965 24 3 5106	sales@trialphaco.com www.trialphaco.com
Latvia	Oriola Riga SIA	Dzelzavas iela 120M Rīga 1021, Latvia	Phone +371 6780 2450 Fax +371 6780 2460	riga@oriola.com www.oriola.lv
Lebanon	Al Yamama	Al Saif Medical Centre Ibn Khaldoun Street Amman, Jordan	Phone +962 6 4655 166 Fax +962 6 4655 167	modhafar@alyamamagroup.com www.alyamamagroup.com
Lithuania	UAB Oriola	Laisvės pr. 75 Vilnius 2022, Lithuania	Phone +370 526 88453 Fax +370 526 88400	giedrius.marcinkonis@oriola.lt www.oriola.com
Luxembourg	Miltenyi Biotec B.V.	Postbus 85183 3508 AD Utrecht The Netherlands	Phone (Lux.) 800 24971 Fax (Lux.) 800 24984	macs@miltenyibiotec.nl www.miltenyibiotec.com

In the United States CliniMACS products for clinical use are available only under an approved Investigational Device Exemption (IDE).

Distributors: products for clinical use

Malaysia	T-Medic Sdn Bhd	H-0-11, Jalan Sri Hartamas 1 (1/70A), 60, Plaza Damas Sri Hartamas, Taman Sri Hartamas 50480 Kuala Lumpur, Malaysia	Phone +60 3 6201 9060 Fax +60 3 6201 9062	www.transmedic.com.sg
Mexico	UNIPARTS, S.A. DE C.V.	Galileo 92, Colonia Polanco 11550 Mexico, D.F., Mexico	Phone +52 55 5281 4718 Fax +52 55 5281 4722	uniparts@uniparts.com.mx www.uniparts.com.mx
Netherlands	Miltenyi Biotec B.V.	Postbus 85183 3508 AD Utrecht The Netherlands	Phone (NL) 0800 4020 120 Fax (NL) 0800 4020 100	macs@miltenyibiotec.nl www.miltenyibiotec.com
New Zealand	Miltenyi Biotec, Australia Pty. Ltd.	Unit 16 A, 2 Eden Park Drive North Ryde, NSW 2113, Australia	Phone +61 02 8877 7400 Fax +61 02 9889 5044	macs@miltenyibiotec.com.au www.miltenyibiotec.com
Northern Ireland	Ulster Anaesthetics Ltd	Maryland Industrial Estate 286 Ballygowan Road Moneyrea BT23 6BL Northern Ireland	Phone +44 2890 448800 Fax +44 2890 449400	vector@ua-ltd.co.uk
Oman	Al Yamama	Al Saif Medical Centre, Ibn Khaldoun Street Amman, Jordan	Phone +962-6-4655 166 Fax +962-6-4655 167	modhafar@alyamamagroup.com www.alyamamagroup.com
Pakistan	T & A Scientific	5, 206-C, 3rd Floor Shahida Mension Central Commercial Area P.E.C.H.S, Block 2 Karachi 75400, Pakistan	Phone +92 21 4555 351 Fax +92 21 4532 574	tascientific@yahoo.com
Peru	Inmunochem	Luis F. Villaran 904 27 Lima, Peru	Phone +51 1 440 9678 Fax +51 1 422 3701	lfquiroz@terra.com.pe www.inmunochem.com
Poland	MEDianus Sp. z o.o.	Ul. Cystersów 13/2 31-553 Kraków, Poland	Phone +48 12 665 3131 Fax +48 12 665 3132	medianus@medianus.net www.medianus.net
Portugal	Citomed Lda.	Rua D. Luis de Noronha no 4, 7º piso 1050-072 Lisboa, Portugal	Phone +351 218 421 430 Fax +351 218 421 439	info@citomed.pt www.citomed.pt
Qatar	Al Yamama	Al Saif Medical Centre, Ibn Khaldoun Street Amman, Jordan	Phone +962 6 4655 166 Fax +962 6 4655 167	modhafar@alyamamagroup.com www.alyamamagroup.com
Romania	NOVAINTERMED	No. 5A, Drumul Potcoavei str. 077190 Pipera-Tunari, Ilfov, Romania	Phone +40 21 230 74 30 Fax +40 21 230 71 30	customer_care_rep2@novaintermed.ro www.novaintermed.ro
Russia	POCARD Ltd.	3-rd Cherepkovskaya 15a 121552 Moscow, Russia	Phone +7 495 414 68 15 Fax +7 495 414 68 20	pocard@cardio.ru www.pocard.ru
Saudi Arabia	Global Health Company	PO. Box 92690 11663 Riyadh, Saudi Arabia	Phone +966 1 4125 252 Fax +966 1 4124 141	feky1@gawab.com
Singapore	Miltenyi Biotec Asia Pacific Pte. Ltd.	100 Beach Road, #28-06 to 28-08 Shaw Tower Singapore 189702	Phone +65 6238 8183 Fax +65 6238 0302	clanimacs@miltenyibiotec.com.sg www.miltenyibiotec.com
Slovak Republic	BIOHEM spol.s.r.o.	Zlátovská 2211 911 01 Trenčín, Slovakia	Phone +421 32 6524 998 Fax +421 32 6525 792	biohem@biohem.sk www.biohem.sk
South Africa	BIOCOM Biotech	PO BOX 13004 0014 Clubview (Centurion) South Africa	Phone +27 82 484 7934 Fax +27 12654 4614	info@biocombiotech.co.za www.biocombiotech.co.za

Distributors: products for clinical use

South Korea	Medilab Korea Co. Ltd.	4 F. WonHyun Building 190-15 Bangyi-Dong, Songpa-Gu Seoul 138-050, Korea	Phone +82 2 424 6367 Fax +82 2 412 6535	clinimacs@medilab.co.kr www.medilabkorea.com
Spain	Miltenyi Biotec S.L.	C/ Luis Buñuel 2 Ciudad de la Imagen 28223 Pozuelo de Alarcón (Madrid), Spain	Phone +34 91 512 1290 Fax +34 91 512 1291	macs@miltenyibiotec.es www.miltenyibiotec.com
Switzerland	Miltenyi Biotec GmbH	Friedrich-Ebert-Straße 68 51429 Bergisch Gladbach, Germany	Phone +49 2204 8306-0 Fax +49 2204 85197	macs@miltenyibiotec.de www.miltenyibiotec.com
Taiwan	Miltenyi Biotec Asia Pacific Pte Ltd.	100 Beach Road, #28-06 to 28-08 Shaw Tower Singapore 189702	Phone +65 6238 8183 Fax +65 6238 0302	clinimacs@miltenyibiotec.com.sg www.miltenyibiotec.com
Thailand	Venus Technology Co., Ltd	636/2 Soi Ramkhamhaeng 39, Theplila 1, Prachautit Road, Wangthonglang, Bangkok 10310, Thailand	Phone +66 2 9356 761 Fax +66 2 9356 766	interco@blood.co.th
Turkey	MAK Sağlık Ürünleri İth. İhr. Tic. San. İtd. Şti.	Cevizlidere Mah. Ceyhun Atuf Kansu cad. No: 137/12 6450 Balgat / Ankara, Turkey	Phone +90 216 360 62 74 Fax +90 216 360 62 75	mak@makmedikal.com www.makmedikal.com
United Arab Emirates	Al Yamama	Al Saif Medical Centre Ibn Khaldoun Street Amman, Jordan	Phone +962 6 4655166 Fax +962 6 4655167	modhafar@alyamamagroup.com www.alyamamagroup.com
Ukraine	MAX-WELL SOCPs	100 Shevchenko str 08300 Boryspil	Phone +38 044 503 60 25 Fax +38 044 503 60 29	d.tatarkin@maxwell-pharma.com www.maxwell-pharma.com
United Kingdom	Miltenyi Biotec Ltd.	Almac House Church Lane, Bisley Surrey GU24 9DR, UK	Phone +44 1483 799 800 Fax +44 1483 799 811	macs@miltenyibiotec.co.uk www.miltenyibiotec.com
United States	Miltenyi Biotec Inc.	12740 Earhart Avenue Auburn, CA 95602, USA	Phone +1 530 888 8871 Fax +1 530 745 2899	macs@miltenyibiotec.com www.miltenyibiotec.com
Venezuela	Biologia Molecular y Biotecnologia	Urbanizacion Terrazas del Club Hipico, Complejo Residencial Brisas del Prado, Terraza A, Centro Comercial Brisas del Prado, Local Comercial No 7, Caracas, Edo Miranda	Phone + 58 212 975 2617 Fax +58 212 978 4180	miguelbio@cantv.net www.biochrom.com.ve
Other countries	Miltenyi Biotec GmbH	Friedrich-Ebert-Straße 68 51429 Bergisch Gladbach, Germany	Phone +49 2204 8306-0 Fax +49 2204 85197	macs@miltenyibiotec.de www.miltenyibiotec.com

CliniMACS® Reagents, Tubing Sets, Instruments and PBS/EDTA Buffer are manufactured and controlled under an ISO 13485 certified quality system. In Europe, the CliniMACS® System and Reagents are available as CE-marked medical devices.

In the USA, CliniMACS® System components including the CliniMACS® Reagents are available for use only under an approved Investigational New Drug (IND) application or Investigational Device Exemption (IDE). CliniMACS® MicroBeads are for research use only and not for use in humans.

Unless otherwise specifically indicated, Miltenyi Biotec products and services are for research only and not for therapeutic or diagnostic use.

MACS and CliniMACS are registered trademarks of Miltenyi Biotec GmbH.

Printed in May 2008

© 2008 Miltenyi Biotec GmbH



Miltenyi Biotec

**Germany/Austria/
Switzerland**

Miltenyi Biotec GmbH
Friedrich-Ebert-Straße 68
51429 Bergisch Gladbach
Germany
Phone +49 2204 8306-0
Fax +49 2204 85197
macs@miltenyibiotec.de

USA/Canada

Miltenyi Biotec Inc.
12740 Earhart Avenue
Auburn, CA 95602, USA
Phone 800 FOR MACS
Phone +1 530 888 8871
Fax +1 530 888 8925
macs@miltenyibiotec.com

Australia

Miltenyi Biotec
Australia Pty. Ltd.
Unit 16A, 2 Eden Park Drive
North Ryde, NSW 2113
Australia
Phone +61 02 8877 7400
Fax +61 02 9889 5044
macs@miltenyibiotec.com.au

Benelux

Miltenyi Biotec B.V.
Postbus 85183, 3508 AD Utrecht
The Netherlands
macs@miltenyibiotec.nl
Customer service Netherlands
Phone 0800 4020120
Fax 0800 4020100
Customer service Belgium
Phone 0800 94016
Fax 0800 99626
Customer service Luxembourg
Phone 800 24971
Fax 800 24984

China

Miltenyi Biotec Trading
(Shanghai) Co., Ltd.
Shanghai Office
Fareast International Plaza A
Rm. 2301, No. 319 Xianxia Rd.
Shanghai 200051, P.R. China
Phone +86 21 62351005
Fax +86 21 62350953
macs@miltenyibiotec.com.cn

France

Miltenyi Biotec SAS
10 rue Mercœur
75011 Paris, France
Phone +33 1 56 98 16 16
Fax +33 1 56 98 16 17
macs@miltenyibiotec.fr

Italy

Miltenyi Biotec S.r.l.
Via Persicetana, 2/D
40012 Calderara di Reno (BO)
Italy
Phone +39 051 6 460 411
Fax +39 051 6 460 499
macs@miltenyibiotec.it

Japan

Miltenyi Biotec K.K.
Nittsu-Eitai Building 5F
16-10 Fuyuki, Koto-ku, Tokyo
135-0041, Japan
Phone +81 3 5646 8910
Fax +81 3 5646 8911
macs@miltenyibiotec.jp

Singapore

Miltenyi Biotec
Asia Pacific Pte Ltd.
100 Beach Road
#28-06 to 28-08 Shaw Tower
Singapore 189702
Phone +65 6238 8183
Fax +65 6238 0302
macs@miltenyibiotec.com.sg

Spain

Miltenyi Biotec S.L.
C/Luis Buñuel 2
Ciudad de la Imagen
28223 Pozuelo de Alarcón
(Madrid), Spain
Phone +34 91 512 12 90
Fax +34 91 512 12 91
macs@miltenyibiotec.es

United Kingdom

Miltenyi Biotec Ltd.
Almac House, Church Lane
Bisley
Surrey GU24 9DR
United Kingdom
Phone +44 1483 799 800
Fax +44 1483 799 811
macs@miltenyibiotec.co.uk