PCR Workstation

PCR1000

Page Finder

Important Information	ii
Waste Electrical & Electronic Equipment (WEEE)	vii
Safety Information	viii
Specifications	ix
Using the PCR Workstation.....	1
Installation and Operation of the PCR Workstation ...	3
Disposal.....	5
Warranty	6
Warning	6
Ordering Information.....	7

Important Information – English

- If this equipment is used in a manner not specified by Hoefer, Inc. the protection provided by the equipment may be impaired.
- This instrument is designed for indoor laboratory use only.
- Only accessories and parts approved or supplied by Hoefer, Inc. may be used for operating, maintaining, and servicing this product.
- Warning! Because this instrument can develop sufficient voltage and current to produce a lethal shock, care must be exercised in its operation.
- This instrument is designed in accordance with the EN61010-1:2001 electrical safety standard. Nevertheless, it should be used only by properly trained operators. Read this entire manual before using the instrument and use only according to the instructions.
- The instrument must always be used with the earth lead of the power cord correctly grounded to earth at the mains outlet.
- Use only undamaged electrical wire and equipment specific for the voltages you will use. All equipment connected to high voltage should be in accordance with EN61010-1:2001.
- Keep the instrument as dry and clean as possible. Wipe regularly with a soft, damp cloth. Let the instrument dry completely before use.
- Do not operate the instrument in extreme humidity (above 80%). Avoid condensation by letting the unit equilibrate to ambient temperature when taking the instrument from a colder to a warmer environment.
- To permit sufficient cooling, ensure that the vents of the instrument are not covered.

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Waste Electrical and Electronic Equipment (WEEE)

English

This symbol indicates that the waste of electrical and electronic equipment must not be disposed as unsorted municipal waste and must be collected separately. Please contact an authorized representative of the manufacturer for information concerning the decommissioning of your equipment.

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Safety Information

These units are capable of delivering potentially lethal voltage when connected to a power supply and are to be operated only by qualified technically trained personnel.

Please read the entire operator's manual thoroughly before operating this unit.

These units comply with the statutory EC safety requirements:

98/79/EC IN VITRO DIAGNOSTIC DIRECTIVE of October 27, 1998

73/23/EEC: LOW VOLTAGE DIRECTIVE:
IEC 1010-1:1990 plus amendment 1:1992

EN 61010-1:1993/BS EN 61010-1:1993

89/336/EC: The EMC directive for electromagnetic compatibility

These units are RoHS compliant.

Hoefer PCR Workstations are designed to give long service and reproducible results in your laboratory. A few moments spent reading these instructions will ensure that your expectations are reflected in the successful use of the apparatus.

First check that the apparatus has been received complete and undamaged following shipment. Any faults or losses must be reported to hoefer immediately. Hoefer cannot accept responsibility for goods returned without prior notification.

Refer to the packing list and check that all components and accessories are present.

Please retain all packaging and materials until the warranty period has expired.

Specifications

Application: For laboratory use only.

Country of Origin: USA

Manufacturer:

Hoefer, Inc.

84 October Hill Road

Holliston, MA 01746

Toll Free: 1-800-227-4750

Phone: 1-508-893-8999

Fax: 1-508-429-5732

E-mail: support@hoeferinc.com

Construction:

Size (h × w × d): 770 × 560 × 420 mm (external)
750 × 540 × 400 mm (internal)

Acrylic Thickness: 10 mm; confers protection
against the UV wavelength and
high energy β -emitters

Mains Supply: 105–250 V AC, 50–60 Hz

Mains Fuse: 2.5 Amp

White Light: 1 × 15 Watt; visible wavelength

UV-C Light: 4 × 15 Watt; 254 nm wavelength

Timer Control: 0–120 minutes

Weight: 19 Kg

Environmental Conditions

- This apparatus is intended for indoor use only.
- This apparatus can be operated safely at an altitude of 2000 m.
- The normal operating temperature range is between 40 °C and 65 °C.
- Maximum relative humidity 80% for temperatures up to 31 °C decreasing linearly to 50% relative humidity at 40 °C.
- The apparatus is rated POLLUTION DEGREE 2 in accordance with IEC 664. POLLUTION DEGREE 2, states that: “Normally only non-conductive pollution occurs. Occasionally, however, a temporary conductivity caused by condensation must be expected”.

UV Safety

The cabinets are manufactured from a material impermeable to UV light. Safety switches on the doors will instantly turn off the UV lights when the doors are opened, however, it is recommended that safety goggles are worn when the UV lights are on as UV-radiation can irreversibly damage eyes and skin.

Using the PCR Workstation

Introduction

The Hoefer[®] PCR Workstation provides effective decontamination of reagents and equipment before carrying out sensitive PCR reactions, particularly when amplifying DNA sequences which are either in limited supply or of low copy number. Four timer-controlled 16-Watt UV bulbs enable the user to control the exposure time and dose of high energy UV irradiation required to denature nucleic acids that cause unwanted background contamination, while a timer override switch allows the UV source to remain on indefinitely if preferred. Safety interlocks switch-off the UV bulbs automatically when the cabinet side doors are opened, preventing accidental exposure to the UV source. The cabinet's construction from UV impermeable 10 mm acrylic also serves as an effective barrier against radioactive isotopes, allowing the user to work in complete safety with high energy β -emitters such as ³²P. A single white light bulb illuminates the work area when the cabinet is in use.

Benefits Include

1. **Four 16-Watt UV bulbs** — generate sufficient high energy UV irradiation to denature nucleic acids in as little as 30 minutes, minimizing unwanted background contamination in PCR reactions.
2. **Optimal reflectors** — provide uniform irradiation of the whole work surface area within the cabinet.
3. **Timer-control** — affords the user total control over exposure time and UV-dosage.
4. **Timer override switch** — for constant UV irradiation.
5. **Safety interlocks** — immediately cut out the UV source when the side doors are opened, preventing accidental UV exposure.
6. **10 mm optical acrylic** — acts as a safety barrier to UV and α and β radiation, allowing the cabinet to be used as a radiation work station if preferred.
7. **White light** — illuminates the work area when the cabinet is in use.
8. **Recommended for use with the RPP-TY6854 safety tray and RPP-TL6854 easy-clean tray liners** (not included).

Installation and Operation of the PCR Workstation

A. Installation

1. Check the cabinet and power cord for any damage. **DO NOT** use if the system shows any sign of damage.
2. Place the cabinet on a flat surface, using a safety tray if required.
3. Connect the power cord into the back of the cabinet and insert into an grounded mains socket.

B. Using the UV Lights

1. Place the objects to be sterilized inside the cabinet and close the side doors.
2. Switch the 'Power' switch to the ON position.
3. Turn the timer to the required time. Twenty minutes is sufficient time to sterilize most pieces of equipment. The UV lights will automatically turn on and then switch off after the set time.
4. If the UV lights are required to remain permanently on, then switch the UV light switch to the ON position. This will override the timer.

C. Using the White Lights

1. Switch the 'Power' switch to the ON position.
2. Switch the 'White Light' switch to the ON position.

Note: The UV lights will not operate when the doors are open.

Note: The white lights are not operated by the timer switch and will remain on when the doors are open.

IMPORTANT: Acrylic plastic is NOT resistant to aromatic or halogenated hydrocarbons, ketones, esters, alcohol's (over 25%) and acids (over 25%). They should NOT be used for cleaning.

CAUTION: The lamps become very hot during use. Make sure that the lamps have cooled down before removing them.

D. General Care and Maintenance

- The unit should be disconnected from the power source prior to cleaning.
- Before use, clean and dry the apparatus with **DISTILLED WATER ONLY.**

DO NOT use abrasive pads, cleaners, creams or scourers. Dry components with clean tissues prior to use. The cabinet should not be used on or above surfaces cleaned with aromatic or halogenated hydrocarbons, ketones, esters, alcohol's (over 25%) and acids (over 25%) or where such compounds have been used because of the risk of vapors damaging the acrylic.

There are no internal operator serviceable parts in this unit. If the power should fail, the unit must be returned to the authorized service center.

E. Replacing the Lamps

Over time, the lamps will diminish in their output. It is therefore recommended that all the lamps are changed at the same time, for uniform efficiency. Please use a qualified safety engineer to replace and check the operation of the lamps.

1. Remove the power cord from the back of the unit.
2. To remove the lamps open both safety doors fully and using both hands, twist the lamps one quarter turn. The lamps will then come out of their sockets.
3. To insert new lamps, place the lamps in their sockets and twist them until they fix in position.
4. Replace the power cord and have the cabinet checked by a qualified safety engineer.

Disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that it must not be disposed of with unsorted waste.

Instead it is your responsibility to dispose of your equipment correctly at lifecycle-end by handing it over to an authorized facility for separate collection and recycling.

It is also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so that the persons involved in the disposal and recycling of the equipment are protected from any potential hazard.

For more information about where you can drop off your waste equipment, please contact your local dealer from whom the equipment was purchased originally.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

Warranty

Hoefer guarantees that the unit you have received has been thoroughly tested and meets its published specification.

This unit (excluding all accessories) is warranted against defective material and workmanship for a period of twelve (12) months from the date of shipment ex factory.

Hoefer will repair all defective equipment returned during the warranty period without charge, provided the equipment has been used under normal laboratory conditions and in accordance with the operating limitations and maintenance procedures outlined in this instruction manual and when not having been subject to accident, alteration, misuse or abuse.

No liability is accepted for loss or damage arising from the incorrect use of this unit. Hoefer's liability is limited to the repair or replacement of the unit or refund of the purchase price, at Hoefer's option. Hoefer is not liable for any consequential damages.

Hoefer reserves the right to alter the specification of its products without prior notice. This will enable us to implement developments as soon as they arise.

Hoefer products are for research use only.

A return authorization must be obtained from Hoefer before returning any product.

Warning

DO NOT attempt to remove the outer casing or make repairs to this or any Hoefer electrical products, should any unit fail.

Contact Hoefer if the need for repair or service is required.

Ordering Information

Product	Qty.	Catalog No.
PCR Workstation UV Sterilization Cabinet for 115 VAC power supply	1	PCR1000-115V
UV Sterilization Cabinet for 230 VAC power supply	1	PCR1000-230V

Accessories

UV Germicidal Bulb, 15 W, 254 nm	1	PCR-UVLS
Radiation Safety Tray, 68 × 54 cm (W × L)	1	RPP-TY6854
Liners, 68 × 54 cm (W × L)	25/pk	RPP-TL6854

Hoefler, Inc.

84 October Hill Road
Holliston, MA 01746

Toll Free: 1-800-227-4750

Phone: 1-508-893-8999

Fax: 1-508-893-0176

E-mail: support@hoeflerinc.com

Web: www.hoeflerinc.com

Hoefler is a registered trademark of
Hoefler, Inc.

© 2012 Hoefler, Inc. —
All rights reserved.

Printed in the USA.