RayBio[®] Human/Mouse/Rat Glucagon Enzyme Immunoassay Kit

Catalog #: EIA-GLU, EIAM-GLU, EIAR-GLU

User Manual Last revised December 1, 2015

Caution: Extraordinarily useful information enclosed



ISO 13485 Certified

3607 Parkway Lane, Suite 100 Norcross, GA 30092

Tel: 1-888-494-8555 (Toll Free) or 770-729-2992, Fax:770-206-2393 Web: www.RayBiotech.com, Email: info@raybiotech.com

Table of Contents

Section				
I.	Introduction			
II.	General Description			
III.	How It Works	4		
IV.	Storage	5		
V.	Reagents			
VI.	Additional Materials Required			
VII.	Reagent Preparation A. Preparation of Plate and Anti-Glucagon Antibody B. Preparation of Biotinylated Peptide (Item F) C. Preparation of Standards D. Preparation of Positive Control E. Preparation of Samples F. Preparation of Wash Buffer and HRP-Strep			
VIII.	Assay Procedure	10		
IX.	Assay Procedure Summary	11		
X.	Calculation of Results A. Typical Data B. Sensitivity C. Detection Range D. Reproducibility E. Assay Diagram	12 12 12 12 12 12 13		
XI.	Specificity			
XII.	Select Publications	14		
XIII.	Troubleshooting Guide			

Please read the entire manual carefully before starting your experiment

I. Introduction

Glucagon is a 29-amino acid peptide hormone secreted by the pancreas. Its function is to raise blood glucose levels to opposing the effect of insulin, which lowers blood glucose levels.

Glucagon is synthesized and secreted from alpha cells of the endocrine portion of the pancreas. In rodents, the alpha cells are located in the outer rim of the islet. However, alpha cells in human pancreas are distributed throughout the islet.

Glucagon and insulin are part of a feedback system that keeps blood glucose levels at a stable level. The pancreas releases glucagon when glucose levels fall too low. Glucagon causes the liver to convert stored glycogen into glucose, which is released into the bloodstream. Glucagon raises blood glucose levels. High blood glucose levels stimulate the release of insulin. Insulin allows glucose to be taken up and used by insulin-dependent tissues.


Glucagon has important clinical applications. Abnormally-elevated levels of glucagon may be caused by pancreatic tumors, such as glucagonoma, with symptoms including necrolytic migratory erythema, reduced amino acids, and hyperglycemia. It may occur alone or in the context of multiple endocrine neoplasia type 1 (MENI).

II. General Description

The RayBio[®] Glucagon Enzyme Immunoassay (EIA) Kit is an in vitro quantitative assay for detecting Glucagon peptide based on the competitive enzyme immunoassay principle.

In this assay, a biotinylated Glucagon peptide is spiked into the samples and standards. The samples and standards are then added to the plate, where the biotinylated Glucagon peptide competes with endogenous (unlabeled) Glucagon for binding to the anti-Glucagon antibody. After a wash step, any bound biotinylated Glucagon then interacts with horseradish peroxidase (HRP)-streptavidin, which catalyzes a color development reaction. The intensity of the colorimetric signal is directly proportional to the amount of captured biotinylated Glucagon peptide and inversely proportional to the amount of endogenous Glucagon in the standard or samples. A standard curve of known concentration of Glucagon peptide can be established and the concentration of Glucagon peptide in the samples can be calculated accordingly.

III. How It Works

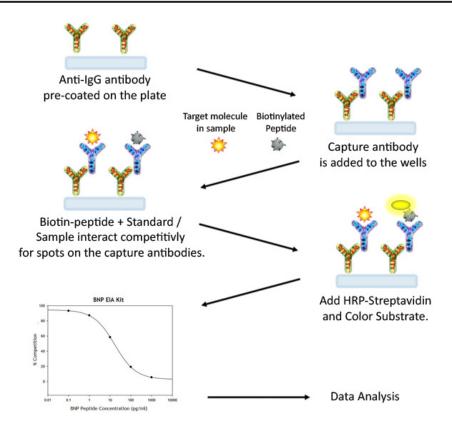

IV. Storage

The entire kit may be stored at -20°C to -80°C for up to 6 months from the date of shipment. For extended storage, it is recommended to store at -80°C. **Avoid repeated freeze-thaw cycles.** For prepared reagent storage, see table below.

V. Reagents

Component	Size / Description	Storage / Stability After Preparation		
Glucagon Microplate (Item A)	96 wells (12 strips x 8 wells) coated with secondary antibody.	1 month at 4°C*		
Wash Buffer Concentrate (20X) (Item B)	25 ml of 20X concentrated solution.	1 month at 4°C		
Standard Glucagon Peptide (Item C)	2 vials of Lyophilized Glucagon Peptide. 1 vial is enough to run each standard in duplicate.	Do not store and reuse		
Anti-Glucagon Polyclonal Antibody (Item N)	2 vials of Lyophilized anti-Glucagon.	Do not store and reuse		
5X Assay Diluent B (Item E)	15 ml of 5X concentrated buffer. Diluent for both standards and samples including serum, plasma, cell culture media or other sample types.	1 month at 4°C		
Biotinylated Glucagon Peptide (Item F)	2 vials of Lyophilized Biotinylated Glucagon Peptide, 1 vial is enough to assay the whole plate.	Do not store and reuse		
HRP-Streptavidin Concentrate (Item G)	600 µl 200X concentrated HRP-conjugated streptavidin.	Do not store and reuse		
Positive Control (Item M)	1 vial of Lyophilized Positive Control.	Do not store and reuse		
TMB One-Step Substrate Reagent (Item H)	12 ml of 3,3,5,5'-tetramethylbenzidine (TMB) in buffer solution.	N/A		
Stop Solution (Item I)	8 ml of 0.2 M sulfuric acid.	N/A		

^{*}Return unused wells to the pouch containing desiccant pack, reseal along entire edge.

VI. Additional Materials Required

- 1. Microplate reader capable of measuring absorbance at 450 nm
- 2. Precision pipettes to deliver 2 µl to 1 ml volumes
- 3. Adjustable 1-25 ml pipettes for reagent preparation
- 4. 100 ml and 1 liter graduated cylinders
- 5. Absorbent paper
- 6. Distilled or deionized water
- 7. SigmaPlot software (or other software which can perform four-parameter logistic regression models)
- 8. Tubes to prepare standard or sample dilutions
- 9. Orbital shaker
- 10. Aluminum foil
- 11. Plastic wrap

VII. Reagent Preparation

Keep kit reagents on ice during reagent preparation steps.


A. Preparation of Plate and Anti-Glucagon Antibody

- 1. Equilibrate plate to room temperature before opening the sealed pouch.
- 2. Label removable 8-well strips as appropriate for your experiment.
- 5X Assay Diluent B (Item E) should be diluted 5-fold with deionized or distilled water.
- Briefly centrifuge the anti-Glucagon antibody vial (Item N) and reconsititute with 55 μl of 1X Assay Diluent B to prepare the antibody concentrate. Pipette up and down to mix gently.
- 5. The antibody concentrate should then be diluted 100-fold with 1X Assay Diluent B. This is your anti-Glucagon antibody working solution, which will be used in step 2 of Assay Procedure (Section VIII).

Note: The following steps may be done during the antibody incubation procedure (step 2 of Assay Procedure)

B. Preparation of Biotinylated Glucagon (Item F)

- 5. Briefly centrifuge the vial of Biotinylated Glucagon (Item F) and reconstitute with 20 µl of ddH₂O before use.
- 6. See the image below for proper preparation of Item F. Transfer the entire contents of the Item F vial into a tube containing 10 ml of 1X Assay Diluent B. This is your Working Stock of Item F. Pipette up and down to mix gently. The final concentration of biotinylated Glucagon will be **80 pg/ml**.
 - a. Second Dilution of Item F for Standards: Add 2 ml of Working Stock Item F to 2 ml of 1X Assay Diluent B. The final concentration of biotinylated Glucagon will be **40 pg/ml**.
 - b. Second Dilution of Item F for Positive Control: Add 100 µl of Working Stock Item F to 100 µl of the prepared Positive Control (Item M). (See section D for Positive Control preparation) The final concentration of biotinylated Glucagon will be **40 pg/ml**.
 - c. Second Dilution of Item F for samples: Add 125 μl of Working Stock Item F to 125 μl of prepared sample (see section E for sample preparation). This is a 2-fold dilution of your sample. The final concentration of biotinylated Glucagon will be 40 pg/ml.

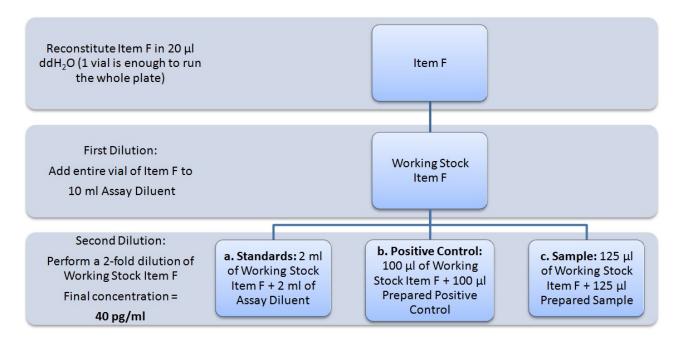


C. Preparation of Standards

7. Label 6 microtubes with the following concentrations: 1000 pg/ml, 100 pg/ml, 10 pg/ml, 0.1 pg/ml and 0 pg/ml. Pipette 450 µl of biotinylated Glucagon Item F working solution (prepapred in step 6a) into each tube, except the 1,000 pg/ml (leave this one empty).

It is very important to make sure the concentration of biotinylated Glucagon is 40 pg/ml in all standards.

- 8. Briefly centrifuge the vial of Glucagon Standard (Item C). Reconstitute with 10 μl of ddH₂O and briefly vortex if desired. Pipette 8 μl of Item C and 792 μl of 40 pg/ml biotinylated Glucagon working solution (prepared in step 6a) into the tube labeled 1000 pg/ml. Mix thoroughly. This solution serves as the first standard (1000 pg/ml Glucagon standard, 40 pg/ml biotinylated Glucagon).
- 9. To make the 100 pg/ml standard, pipette 50 µl of the 1000 pg/ml Glucagon standard into the tube labeled 100 pg/ml. Mix thoroughly.
- 10. Repeat this step with each successive concentration, preparing a dilution series as shown in the illustration below. Each time, use 450 μl of biotinylated Glucagon and 50 μl of the prior concentration until the 0.1 pg/ml is reached. Mix each tube thoroughly before the next transfer.

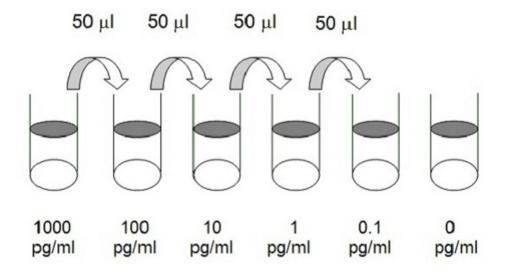

D. Positive Control Preparation

- 11. Briefly centrifuge the Positive Control vial (Item M) and reconstitute with 100 μl of ddH₂O.
- 12. Refer to step 6b. This is a 2-fold dilution of the Positive Control. The final concentration of biotinylated Glucagon should still be 40 pg/ml.

The Positive Control is a cell culture media sample that serves as a system control to verify that the kit components are working. The resulting OD will not be used in any calculations; if no positive competition is observed please contact RayBiotech Technical Support. The Positive Control may be diluted further if desired, but be sure the final concentration of biotinylated Glucagon is 40 pg/ml.

E. Sample Preparation

13. If you wish to perform a 2-fold dilution of your sample, proceed to step 6c. If you wish to perform a higher dilution of your sample, dilute your sample with 1X Assay Diluent B before performing step 6c.

EXAMPLE (to make a 4-fold dilution of sample):

- a. Dilute sample 2-fold (62.5 µl of sample + 62.5 µl of 1X Assay Diluent B.).
- b. Perform step 6c (125 μ l of working solution Item F + 125 μ l of sample prepared above).

The total volume is 250 µl, enough for duplicate wells on the microplate.

It is very important to make sure the final concentration of the biotinylated Glucagon is **40 pg/ml**.

Note: Optimal sample dilution factors should be determined empirically, however you may reference below for recommended dilution factors for serum: Human=2X Mouse=2X Rat=2X.

If you have any questions regarding the recommended dilutions you may contact technical support at 888-494-8555 or techsupport@raybiotech.com.

F. Preparation of Wash Buffer and HRP

- 14. If Item B (20X Wash Concentrate) contains visible crystals, warm to room temperature and mix gently until dissolved.
- 15. Dilute 20 ml of Wash Buffer Concentrate into deionized or distilled water to yield 400 ml of 1X Wash Buffer.
- 16. Briefly centrifuge the HRP-Streptavidin vial (Item G) before use.
- 17. Dilute the HRP-Streptavidin concentrate 200-fold with 1X Assay Diluent B.

VIII. Assay Procedure

- 1. Keep kit reagents on ice during reagent preparation steps. It is recommended that all standards and samples be run at least in duplicate.
- 2. Add 100 μl of Anti-Glucagon Antibody (Item N) (See Reagent Preparation step 3) to each well. Incubate for 1.5 hours at room temperature with gentle shaking (1-2 cycle/sec). You may also incubate overnight at 4°C.
- 3. Discard the solution and wash wells 4 times with 1X Wash Solution Buffer (200-300 µl each). Washing may be done with a multichannel pipette or an automated plate washer. Complete removal of liquid at each step is essential to good assay performance. After the last wash, remove any remaining Wash Buffer by aspirating or decanting. Invert the plate and blot it against clean paper towels.
- 4. Add 100 μl of each standard (see Reagent Preparation Section C), Positive Control (see Reagent Preparation Section D) and sample (see Reagent Preparation Section E) in appropriate wells. Be sure to include a blank well (Assay Diluent only). Cover wells and incubate for 2.5 hours at room temperature with gentle shaking (1-2 cycles/sec) overnight or at 4°C.
- 5. Discard the solution and wash 4 times as directed in Step 3.

- 6. Add 100 μl of prepared HRP-Streptavidin solution (see Reagent Preparation step 7) to each well. Incubate for 45 minutes at room temperature with gentle shaking. It is recommended that incubation time should not be shorter or longer than 45 minutes.
- 7. Discard the solution and wash 4 times as directed in Step 3.
- Add 100 μl of TMB One-Step Substrate Reagent (Item H) to each well. Incubate for 30 minutes at room temperature in the dark with gentle shaking (1-2 cycles/sec).
- 9. Add 50 µl of Stop Solution (Item I) to each well. Read at 450 nm immediately.

IX. Assay Procedure Summary

- 1. Prepare all reagents, samples and standards as instructed.
- 2. Add 100 μl anti-Glucagon to each well. Incubate 1.5 hours at room temperature or overnight at 4°C.
- 3. Add 100 µl standard or sample to each well. Incubate 2.5 hours at room temperature or overnight at 4°C.
- Add 100 μl prepared Streptavidin solution. Incubate 45 minutes at room temperature.
- 5. Add 100 µl TMB One-Step Substrate Reagent to each well. Incubate 30 minutes at room temperature.
- Add 50 μl Stop Solution to each well. Read at 450 nm immediately.


X. Calculation of Results

Calculate the mean absorbance for each set of duplicate stands, controls, and samples and subtract the blank optical density. Plot the standard curve using SigmaPlot software (or other software which can perform four-parameter logistic regression models), with standard concentration on the x-axis and percentage of absorbance (see calculation below) on the y-axis. Draw the best-fit curve through the standard points.

Percentage absorbance = (B-blank OD)/ B_0 -blank OD) where B = OD of sample or standard and B_0 = OD of zero standard (total binding)

A. Typical Data

These standard curves are for demonstration only. A standard curve must be run with each assay.

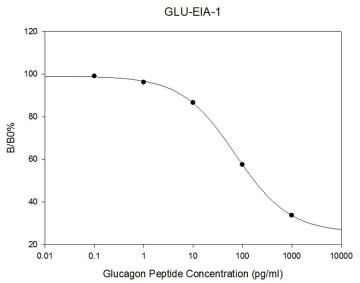

B. Sensitivity

The minimum detectable concentrations of Glucagon is 4.77 pg/ml.

C. Detection Range

1-1,000 pg/ml

D. Reproducibility

Intra-Assay: CV<10% Inter-Assay: CV<15%

E. Assay Diagram

Recommended Plate Layout:

Blank	Blank	SA1	SA1	SA9	SA9	SA17	SA17	SA25	SA25	SA33	SA33
Total Binding	Total Binding	SA2	SA2	SA10	SA10	SA18	SA18	SA26	SA26	SA34	SA34
Standard1	Standard1	SA3	SA3	SA11	SA11	SA19	SA19	SA27	SA27	SA35	SA35
Standard2	Standard2	SA4	SA4	SA12	SA12	SA20	SA20	SA28	SA28	SA36	SA36
Standard3	Standard3	SA5	SA5	SA13	SA13	SA21	SA21	SA29	SA29	SA37	SA37
Standard4	Standard4	SA6	SA6	SA14	SA14	SA22	SA22	SA30	SA30	SA38	SA38
Standard5	Standard5	SA7	SA7	SA15	SA15	SA23	SA23	SA31	SA31	SA39	SA39
Pos Control	Pos Control	SA8	SA8	SA16	SA16	SA24	SA24	SA32	SA32	SA40	SA40

Key:

Blank = Buffer Only

Total Binding = Biotin-Glucagon only

Standard 1 = 1000 pg/ml

Standard 2 = 100 pg/ml

Standard 3 = 10 pg/ml

Standard 4 = 1 pg/ml

Standard 5 = 0.1 pg/ml

Pos Control = Biotin with Item M

XI. Specificity

Cross Reactivity: This EIA kit shows no cross-reactivity with any of the cytokines tested: Ghrelin, Nesfatin, Angiotensin II, NPY and APC

XIV. Publications Citing This Product

 Elliot A., Ustione A., Piston D. Somatostatin and insulin mediate glucose-inhibited glucagon secretion in the pancreatic ?-cell by lowering cAMP. American Journal of Physiology - Endocrinology and Metabolism, 15 January 2015 Vol. 308 no. 2, E130-E143DOI: 10.1152/ajpendo.00344.2014

Species: Mouse

Sample Type: Conditioned Media

XIII. Troubleshooting Guide

Problem	Cause	Solution				
Poor standard curve	Inaccurate pipettingImproper standard dilution	Check pipettes Briefly centrifuge Item C and dissolve the powder thoroughly by gently mixing				
Low signal	 Improper preparation of standard and/or biotinylated antibody Too brief incubation times Inadequate reagent volumes or improper dilution 	 Briefly spin down vials before opening. Dissolve the powder thoroughly. Ensure sufficient incubation time; assay procedure step 2 may be done overnight Check pipettes and ensure correct preparation 				
Large CV	Inaccurate pipettingAir bubbles in wells	Check pipettes Remove bubbles in wells				
High background	 Plate is insufficiently washed Contaminated wash buffer 	 Review the manual for proper wash. If using a plate washer, ensure that all ports are unobstructed. Make fresh wash buffer 				
Low sensitivity	Improper storage of the ELISA kit Stop solution	 Follow storage recomendations in sections IV and V. Keep substrate solution protected from light. Add stop solution to each well before reading plate 				

RayBio[®] ELISA Kits

Over 2,000 ELISA kits available, visit www.RayBiotech.com/ELISA-Kits.html for details.

This product is for research use only.



©2015 RayBiotech, Inc