

# UVL-4F Longwave UV Lamp

## User's Manual

**Important: Please read the instructions carefully prior to using the lamp in order to avoid malfunction or damage.**

### Operation (Fig. 1)

**Spotlight:** Switch the lamp to the | position.


**Fluorescent Lamp:** Turn the fluorescent knob located on the top counter-clockwise.

### Replacing the Fluorescent Tube (Fig. 2)


1. Remove the batteries.
2. Turn the fluorescent knob counter-clockwise until the fluorescent lamp cover is fully exposed.
3. Lift the fluorescent lamp cover up and pull it out.
4. Turn the tube in either direction and replace with a new 6" four-watt fluorescent tube (see replacement parts for ordering information).
5. Put the lamp cover back to the original position.

### Replacing the Bulb (Fig. 2)

1. Turn the spotlight cover counter-clockwise and remove it.
2. Loosen the bulb by using the bulb remover key (key shown in Fig. 1) to turn it counter-clockwise.
3. Put the reflector and the spotlight cover back to the original position.

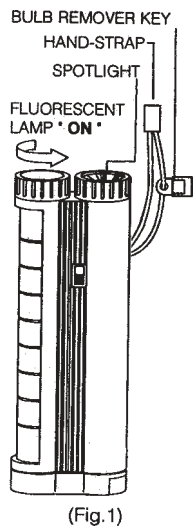

(Fig.1)

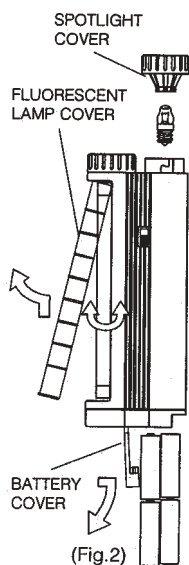

(Fig.2)

### Replacing the Batteries (Fig. 2)

1. Slide the battery cover out and lift it up.
2. Remove the old batteries and replace with four new "AA" size "UM3" batteries, noting the positive and negative indication mark for proper insertion.
3. Slide back the battery cover.

### Maintenance

Keep the lamp away from extreme temperature changes, excessive humidity, moisture, rain or other harsh environments. Do not mix old and new batteries together. It is not recommended to use the fluorescent lamp function when the energy level is low, it will blacken the fluorescent tube. Remove the batteries if the lantern is not to be used for a long period.

### Replacement Parts

	Part Number
Replacement Tube, 4W longwave BLB	34-0010-01



**Corporate Headquarters:** UVP, Inc. 2066 W. 11th Street, Upland, CA 91786 Tel: (800)452-6788 or (909)946-3197 Fax: (909)946-3597 E-Mail: [uvp@uvp.com](mailto:uvp@uvp.com)

**European Sales Operations: Ultra-Violet Products Ltd.** Unit 1, Trinity Hall Farm Estate, Nuffield Road, Cambridge CB4 1TG UK Tel: +44(0)1223-420022 Fax: +44(0)1223-420561 E-Mail: [uvp@uvp.co.uk](mailto:uvp@uvp.co.uk)

