

EyeLink[®] Data Viewer User's Manual

Document Version 2.1.1

Please report all functionality comments and bugs to:

support@sr-research.com

Copyright ©2002-2014, SR Research Ltd.

EyeLink is a registered trademark of SR Research Ltd., Mississauga, Canada

Table of Contents

1	Introduction.....	1
1.1	Data Visualization.....	1
1.2	Event Selection	2
1.3	Interest Area Definition	2
1.4	Event Filtering	2
1.5	Interest Periods and Reaction Time Definition.....	3
1.6	Data Output and Analysis	3
1.7	Experiment Integration	4
2	Installation.....	5
2.1	System Requirements.....	5
2.2	Setup	5
2.2.1	For Standard Installation (applicable to most users).....	5
2.2.2	For Installation using Network Licensing.....	6
3	Working with Files	8
3.1	Creating a Data Viewing Session (DVS).....	8
3.2	Saving a Viewing Session.....	8
3.3	Opening an Existing Viewing Session.....	8
3.4	Saving an Existing Viewing Session to a Different File.....	9
3.5	Importing EyeLink Data Files	9
4	Data Viewer Windows.....	11
4.1	Inspector Window	11
4.1.1	Data Tab.....	11
4.1.1.1	Data Tree Hierarchy.....	13
4.1.1.2	Trial Grouping	13
4.1.1.3	Trial Summary	14
4.1.1.4	Interest Area Templates	16
4.1.1.5	Custom Interest Area Set/Empty Interest Area Set.....	16
4.1.2	Preferences Tab.....	16
4.2	Trial View Window	16
4.2.1	Selecting Data Elements to Display.....	17
4.2.2	Cloning a Trial View	17
4.2.3	Tiling/Cascading Windows.....	17
4.3	Spatial Overlay Trial View	18
4.4	Temporal Graph Trial View.....	20
4.4.1	Viewing Sample and Target Data.....	20
4.4.2	Zooming Operations	22
4.5	Animation View.....	22
5	Working with Events, Samples, and Interest Areas.....	25
5.1	Common Operations	25
5.1.1	Selecting Events.....	26
5.1.2	Deleting Events.....	26
5.2	Fixations.....	26
5.2.1	Selecting Fixations.....	29
5.2.2	Merging Fixations.....	29
5.2.3	Drift Correcting Fixations.....	29

5.2.3.1	Drift-correcting a Group of Selected Fixations	29
5.2.3.2	Precise Drift Correction of a Single Fixation	30
5.2.3.3	Cleaning Fixations	31
5.3	Saccades	32
5.3.1	Merging Saccades	33
5.4	Blink Events	33
5.5	Messages	34
5.5.1	Import Message List	36
5.6	Button Events	37
5.7	Samples	39
5.8	Interest Areas	39
5.8.1	Format of static and dynamic interest areas	40
5.8.2	Interest Area Properties	43
5.8.3	Manually Creating Interest Areas	45
5.8.4	Creating Interest Area: Image Segmentation	48
5.8.5	Saving Interest Areas	48
5.8.6	Importing Interest Area File (for a specific trial)	48
5.8.7	Emptying an Interest Area Set	49
5.8.8	Loading Interest Area Templates	49
5.8.9	Using Interest Area Templates	50
5.8.10	Deleting Interest Area Templates	51
5.8.11	Selecting Interest Areas	51
5.8.12	Copying and Pasting Interest Areas	51
5.8.13	Cleaning Up Dynamic Interest Areas	52
5.8.14	Mouse Record Mode	52
5.8.15	Moving an Interest Area	53
5.8.16	Resizing Rectangular and Elliptical Interest Areas	54
5.8.17	Resizing Freehand Interest Areas	54
5.8.18	Adding a Point to a Freehand Interest Area	54
5.8.19	Deleting an Existing Point from a Freehand Interest Area	55
5.8.20	Deleting Interest Areas	55
5.8.21	Splitting Interest Areas	55
5.8.22	Merging Interest Areas / Instances of a Dynamic Interest Area	56
5.8.23	Conversion between a Static Interest Area and a Dynamic Interest Area ..	56
5.8.24	Examples of Creating Dynamic Interest Areas in a Viewing Session	56
5.8.25	Output Reports and Preferences	58
6	Data Analysis and Output	60
6.1	Managing Trial Variables	60
6.1.1	Trial Variable Manager	61
6.1.2	Trial Variable Value Editor	61
6.1.3	Trial Grouping	62
6.2	Fixation Map	63
6.3	Managing Reaction Time Definitions	68
6.3.1	Reaction Time Manager	69
6.3.2	Viewing Reaction Time Information	72
6.4	Interest Periods	73

6.4.1	Creating Interest Periods.....	73
6.4.2	Using Interest Period.....	75
6.5	Trial Report.....	76
6.5.1	Obtaining a Trial Report.....	76
6.5.2	Variables in the Trial Report.....	78
6.5.3	Using Trial Report.....	81
6.5.4	Configuring the Trial Report.....	82
6.6	Fixation Report.....	83
6.6.1	Obtaining a Fixation Report.....	83
6.6.2	Variables in the Fixation Report.....	85
6.6.3	Configuring the Fixation Report.....	91
6.7	Saccade Report.....	92
6.7.1	Obtaining a Saccade Report.....	92
6.7.2	Variables in the Saccade Report.....	92
6.8	Interest Area Report.....	93
6.8.1	Obtaining an Interest Area Report.....	93
6.8.2	Variables in the Interest Area Report.....	95
6.8.3	Using Interest Area Report.....	101
6.8.3.1	Fixation Sequence Analysis.....	102
6.8.3.2	Regression Analysis.....	103
6.9	Sample Output Report.....	105
6.9.1	Loading Samples into a Viewing Session.....	105
6.9.2	Obtaining a Sample Output Report.....	105
6.9.3	Variables in the Sample Output Report.....	107
6.10	Message Output Report.....	111
6.10.1	Obtaining a Message Output Report.....	112
6.10.2	Variables in the Message Output Report.....	113
6.11	Recording Event Sequence Data.....	116
6.11.1	Obtaining a Recording Event Sequence Data.....	116
6.11.2	Formats of the Recording Event Sequence Data.....	116
6.11.2.1	Fixation Events.....	117
6.11.2.2	Saccade Events.....	118
6.11.2.3	Blink Events.....	118
6.11.2.4	Message Events.....	118
6.11.2.5	Button Events.....	118
7	Protocol for EyeLink Data to Viewer Integration.....	119
7.1	Defining the Start and End of a Trial.....	119
7.1.1	Trial Start Message.....	119
7.1.2	Trial End Message.....	120
7.2	Pre-Trial Message Commands.....	120
7.2.1	Trial Variable Labels.....	120
7.2.2	Display Coordinates.....	121
7.2.3	Trial Grouping.....	121
7.3	Trial Message Commands.....	122
7.3.1	Trial Variable Values.....	122
7.3.2	Single Trial Variable Message Token.....	122

7.3.3	Image Commands	123
7.3.3.1	Image Loading – Fill Full Screen	123
7.3.3.2	Image Loading – Top Left	124
7.3.3.3	Image Loading – Centered	124
7.3.4	Simple Drawing	125
7.3.4.1	Clear Overlay View	125
7.3.4.2	Line drawing	126
7.3.4.3	Drawing a Rectangle	126
7.3.4.4	Drawing a Filled Rectangle	127
7.3.4.5	Drawing a Fixation Point	127
7.3.4.6	Draw List File	128
7.3.5	Interest Area Commands	129
7.3.5.1	Rectangular Interest Area	129
7.3.5.2	Elliptical Interest Area	129
7.3.5.3	Freehand Interest Area	130
7.3.5.4	Interest Area Set	130
7.3.6	Target Position Commands	131
7.4	Reaction Time Definitions	132
7.4.1	Button Events	132
7.4.2	Fixation Events:	133
7.4.3	Saccade Events:	134
7.4.4	Message Events:	135
7.4.5	Reaction Time Definition Set	135
7.5	Changes in the Sample Experiments	136
7.5.1	“Simple” Template	136
7.5.2	“Text” Template	136
7.5.3	“Picture” Template	137
7.5.4	“EyeData” Template	137
7.5.5	“GCWindow” Template	138
7.5.6	“Control” Template	138
7.5.7	“Dynamic” Template	139
7.5.8	“Comm_simple” Template	139
7.5.9	Other Templates	139
8	Preference Settings	140
8.1	General Preferences	141
8.2	Data Loading Preferences	142
8.3	Output / Analysis	144
8.4	Fixation Map	147
8.5	Data Filter Preferences	149
8.6	General Trial View Preferences	152
8.7	Spatial Overlay View Preferences	154
8.8	Image Segmentation Preferences	156
8.9	Temporal Graph Preferences	157
8.10	Animation Preferences	159
9	Release History	160

List of Figures

Figure 4-1. Viewer Display after Loading the Sample Picture.edf File	11
Figure 4-2. Components of the Inspector Window (Left: Data Tab; Right: Preference Tab).....	13
Figure 4-3. Tiling Trial View Windows.	18
Figure 4-4. Sample Spatial Overlay Trial View	19
Figure 4-5. Sample Temporal Graph Trial View.....	20
Figure 4-6. Sample Playback Animation View.	23
Figure 5-1. Sample Spatial Overlay View of Fixation Events.....	27
Figure 5-2. Diagram for Angle Calculation.....	28
Figure 5-3. Drift Correcting Fixations on One Line of Text.	30
Figure 5-4. Options for Fixation Cleaning.....	31
Figure 5-5. Sample Spatial Overlay View of Saccade Events.....	32
Figure 5-6. Sample Temporal Graph View of Blink Events	34
Figure 5-7. Sample Temporal Graph View of Message Events (Note that the default size of the message symbol in the temporal-graph preference settings has been changed).	36
Figure 5-8. Sample Temporal Graph View Showing the Button Events (Note that the default size of the button symbol in the temporal-graph preference settings has been changed).....	38
Figure 5-9. Sample Spatial Overlay View (left) and Temporal Graph View (right) of Raw Sample Data.....	39
Figure 5-10. Sample Text Display Overlaid with Interest Areas.....	40
Figure 5-11. Sample Interest Area Instances in One Interest Period.....	42
Figure 5-12. Sample Interest Area Instances in Another Interest Period	43
Figure 5-13. Creating a Freehand Interest Area.....	46
Figure 5-14. Sample Interest Area Instances in Another Interest Period	47
Figure 5-15. Interest Area Templates Listed in the Tree View	50
Figure 5-16. Mouse Record Mode.....	53
Figure 6-1. Trial Variable Manager	61
Figure 6-2. Trial Variable Value Editor Dialog.....	62
Figure 6-3. Applying Trial Grouping by a Trial Condition Variable and Setting Interest Area Template for a Group of Trials.	63
Figure 6-4. Example fixation map	64
Figure 6-5. Creating a Fixation Map.....	65
Figure 6-6. Types of Fixation Map.....	67
Figure 6-7. Adjusting the Scale of the Fixation Map.....	68
Figure 6-8. Reaction Time Definition Manager	69
Figure 6-9. Reaction Time Definition Editor (General Tab).....	70
Figure 6-10. Reaction Time Definition Editor (Reaction Time Tab)	70
Figure 6-11. Deleting Reaction Time Definitions	72
Figure 6-12. Reaction Time Graphics in Both the Spatial Overlay View and the Temporal Graph View of the Trial View Window.....	72
Figure 6-13. Interest Period Editor (General Tab).....	74
Figure 6-14. Interest Period Editor (Details Tab).....	74
Figure 6-15. Creating a Trial Report.....	76

Figure 6-16. Variables used in the Trial Report.	77
Figure 6-17. Saving the Output Report.	78
Figure 6-18. Using Interest Period for Event Filtering.	82
Figure 6-19. Creating a Fixation Report.	83
Figure 6-20. Variables in a Fixation Report.	84
Figure 6-21. Creating a Saccade Report.	92
Figure 6-22. Variables in a Saccade Report.	93
Figure 6-23. Creating an Interest Area Report.	94
Figure 6-24. Variables in an Interest Area Report.	94
Figure 6-25. Progress Dialog Box Displayed when Creating the Interest Area Report	102
Figure 6-26. Creating a Sample Report.	106
Figure 6-27. Variables in a Sample Report.	106
Figure 6-28. Creating a Message Report.	112
Figure 6-29. Variables in a Message Report.	113
Figure 8-1. Saving properties as defaults.	140

1 Introduction

The EyeLink Data Viewer is a tool that allows the display, filtering, and report output of EyeLink I , EyeLink II, EyeLink 1000, and EyeLink 1000 Plus EDF data files. Data Viewer includes the following areas of functionality.

1.1 Data Visualization

The EyeLink Data Viewer supports three trial-based viewing modes: the Spatial Overlay View, the Temporal Graph View, and the Animation View. In the first two viewing modes, the user can specify which event types to display, including fixations, saccades, blinks, messages, and buttons. Sample traces can also be displayed. The animation viewing mode plays back the gaze data of a selected trial.

The Spatial Overlay View allows a trial to be viewed superimposed on the background image that the participant was looking at while the data was being recorded. This view is ideal for relatively static trial presentations, like reading or visual search paradigms. Fixations can be displayed as either circles or as a scan path. Saccades can be displayed as lines with an arrow specifying the saccade direction. Blinks can be displayed as a line joining the last valid sample position prior to the blink and the first valid sample following the blink. Message and button events can be displayed as a small rectangle, where the position of the event is determined by the position of the nearest sample recorded at the message or button time.

The Temporal Graph View provides a two dimensional plot of trial data. The X axis represents time, while the Y axis represents the location of the samples or events being plotted. This view is suitable for a wide range of experimental paradigms, including those using dynamic trial displays. The scale of the graph can be changed via zoom in and zoom out operations. This view supports binocular display of sample position and pupil size traces.

The color of individual events, as well as the global color of a given event type or sample trace, can be changed as desired. There is a vast number of preferences that can be configured by the user, saved, and then applied to future loaded data.

The Animation View plays back the subject's gaze data in a trial, with a concurrent time code displaying the time from the start of trial recording. If the image that was shown during the trial recording is available, it can be loaded as the background for the playback view. A trackbar control displays the progress of the playback and lets the user select a specific position to start playback. The speed of playback can also be adjusted.

Multiple viewing windows can be opened at one time, displaying data from different trials or displaying data of one single trial in the Spatial Overlay View, Temporal Graph View, and Animation View concurrently.

1.2 Event Selection

EyeLink events can be selected for either of the data views (temporal graph view or spatial overlay view). Detailed properties of the selected event can be examined within a Inspector Window that also provides a list view of all visible events in the selected trial. Properties of a selected event that are editable by the user (like the events color) can be modified directly within the inspector window.

1.3 Interest Area Definition

The Spatial Overlay View allows creation of any number of interest areas for the trial. Rectangular, elliptical, and freehand interest area shapes can be created. The position and shape of the Interest Area can be modified after Interest Area creation.

Users can view fixation-based statistics for a selected interest area, including the fixation count and total dwell time for the interest area, as well as the proportion of fixations and dwell time relative to the trial totals.

Interest area templates can be created which can then be applied to multiple trials, speeding the process of defining interest areas for multiple trials.

1.4 Event Filtering

The EyeLink Data Viewer supports data filtering, including deleting, merging, and drift correcting event types. Not all filtering functions are available to all event types. The original EyeLink EDF file is never modified by the Data Viewer, so the originally recorded data is always available.

Any event can be hidden within the Data Viewer either manually or through the definition of some spatial or temporal filtering preferences. Hiding events allows the user to focus on a subset of a trial's data. Hidden events can be made visible again at a later date if required.

Events can also be deleted from within the Data Viewer. This is useful, for example, if the researcher inspects a specific fixation or saccade event and determines that the event detection performed by the EyeLink on-line parser did not satisfy their needs. If a fixation or saccade event is deleted, the surrounding saccade or fixation events are automatically merged to keep the event structure consistent.

The position of a fixation event can also be adjusted if required. This is only suggested if it is obvious to the user that improper system setup or calibration has resulted in a trial's fixation data containing significant drift that can be easily corrected manually. Any event that is adjusted in this fashion is flagged so that it is clear what events have been manually altered. Optionally, if the position of a fixation is adjusted, the adjoining

saccade end point and start point are also adjusted so that the saccade positions are kept consistent with fixation positions.

1.5 Interest Periods and Reaction Time Definition

For each trial, the user may selectively view and perform analysis on data within a specific period of time within a trial (Interest Period). The EyeLink Data Viewer allows the user to create interest periods based on messages and button events. The created interest periods are added to the interest period list. The user can navigate between different interest periods by selecting the desired interest period from the list. For each interest period, only those events falling within that period are shown in the spatial overlay, temporal graph views, and in the event list of the inspector window. In addition, only those events and data within the selected interest period (custom-defined interest period, or the default – “Full Trial Period”) are exported to the output file.

One important variant of the interest period is the reaction-time period. Reaction Time Definitions (RTD) can be created and applied to all trials loaded into an EyeLink Data Viewer session. Each RTD can have a set of trial condition variables that must be matched for the RTD to be applied to a given trial. The message event to use as the start time for the reaction time can be defined. Each RTD can have a fixation, saccade, button, or message event as the ending event for the reaction time calculation. Depending on the event type selected for the reaction time end event, a set of event properties can be specified to determine which event should be used within each trial. A number of RTDs can be created for a given Data Viewer session, allowing multiple reaction time conditions to be specified. Each trial will only be matched to at most one RTD.

1.6 Data Output and Analysis

Both Spatial Overlay and Temporal Graph trial views can be saved as JPEG images for illustrative purposes.

A series of output reports can be run for a Data Viewer session.

The Fixation Output Report generates one row for each visible fixation within the Data Viewer. A set of output variables can be selected, with each variable output as a column in the report. Fixation report variables include start and end time, duration, average position, interest area identifier, etc. A set of relative variables is also available, including the previous and next fixations position, angle, distance, and direction.

The Saccade Output Report generates one row for each visible saccade within the Data Viewer. A set of output variables can be selected, with each variable output as a column in the report. Saccade output report variables include start and end time, start and end position, amplitude, angle, direction, average and peak velocity, etc. A set of relative variables is also available, including the previous and next saccades start and end position, etc.

The Interest Area Output Report generates one row for each interest area in every trial loaded in the Data Viewer. Variables include the interest area identifier, label, fixation index list, number fixations, summed dwell time, and proportions of fixations and dwell time in the interest area relative to the trial.

The Trial Output Report generates one row of data for each trial within the Data Viewer. Variables include the count of each event type visible within the trial, the average fixation duration and saccadic amplitude, as well as the reaction time calculated for the trial based on the reaction time definitions created for the Data Viewer session.

The Sample Output Report provides a columnar output of eye movement data for each eye sample in a trial. Variables include the index of the sample in a trial, time of the sample, position (x, y), velocity, acceleration, and pupil data of the current sample. If appropriate target position messages are read from the EDF file, the position, velocity, and acceleration data of the targets (up to two targets) can also be provided.

The user can also save the spatial overlay and temporal graph views into image files, save the playback of the trial into a movie file, or create a "landscape" view (i.e., fixation map) for a trial or for a group of trials with the same background image.

1.7 Experiment Integration

The EyeLink Data Viewer allows commands to be inserted directly into the EyeLink EDF file when an experiment is running. This allows the experiment to be significantly integrated into the Data Viewer. These commands are interpreted by the Data Viewer when the data file is loaded and automates several areas of the viewer functionality.

By using a set of predefined viewer commands, the programmer of an experiment can inform the Data Viewer regarding:

- a) The image(s) to load for each trial's Spatial Overlay View.
- b) The Interest Areas to load for each trial.
- c) The Reaction Time Definitions to apply to the Data Viewing session.
- d) The trial condition variables and values to use for each trial.
- e) The position of the target in a dynamic display.

If you use SR Research Experiment Builder software, please make sure you use the Experiment Builder Project Checklist to check up the experiment project before data collection. This checklist is contained in the Experiment Builder User Manual (or from this link <https://www.sr-support.com/forums/showthread.php?t=1012>).

2 Installation

2.1 System Requirements

Operating System: Windows XP, Vista, Windows 7, or Windows 8

Processor: Pentium 4 1.4 GHz or higher (avoid computers with atom CPU)

RAM: Minimum 1GB MB RAM (2GB or higher recommended)

Disk Space: 240 MB free disk space for application plus space required for EyeLink data files and saved viewing sessions.

Graphics Card: video card with Opengl 2.0 support and with at least 128 MB RAM

Monitor Resolution: 1024x768 or better resolution

Other: Free USB port required

2.2 Setup

The latest version of EyeLink Data Viewer installer can be downloaded from <https://www.sr-support.com/forums/showthread.php?t=10>. If you have a previous version of EyeLink Data Viewer installed on the computer, please uninstall it with the DV installer or with Windows® Control Panel tool "Add or Remove Programs" before installing the new version. By default, the EyeLink Data Viewer software will be installed at "{Windows Drive}:\Program Files\SR Research\DataViewer".

Important: A USB Dongle provided by SR Research is required to run this application and must remain connected to the PC while the application is running. If a USB Dongle is not detected by the program, the application will run in a DEMO mode supporting limited data loading.

2.2.1 For Standard Installation (applicable to most users)

The following installation instruction is applicable to users who use a standalone USB dongle that supports a single-PC license.

1. Install the EyeLink Data Viewer software. Double click on the EyeLinkDV1.*.exe installer, keeping the default settings.

2. Install standalone HASP key driver (if this is the first time that the USB dongle have been used on the display PC). You may install the driver by clicking "Start -> All Programs -> SR Research -> Install HASP Driver" from your computer desktop or double clicking on "hdd32.exe" in "C:\Program Files\SR Research\Common" folder.
3. Plug the dongle to the display computer. Click on the License Manager utility (Start -> All Programs -> SR Research -> License Manager) to check for the licensing status for the Data Viewer software.
4. From the computer desktop, click "Start -> All Programs -> SR Research -> EyeLink -> Data Viewer -> DataViewer" to run the application.

2.2.2 For Installation using Network Licensing

The following is applicable to the users who have purchased a network license (i.e., a shared license for several computers on a network that running EyeLink Data Viewer at the same time) for the Data Viewer software.

1. Install the Data Viewer software. Double click on the EyeLinkDV1.*.exe installer, keeping the default settings on InstallShield Wizard screens except for the following two:
 - On the "Setup Type" dialog box, select "Custom".
 - On the "Select Feature" screen, make sure that both "HASP4" and "HASPHL" driver options are selected.
2. Install network HASP key driver. You may install the driver by clicking "Start -> All Programs -> SR Research -> HASP HL Driver" from your computer desktop or double clicking on "HaspUserSetup.exe" in "C:\Program Files\SR Research\Common" folder.
3. Install network HASP License Manager. You may install the tool by clicking "Start -> All Programs -> SR Research -> Networked HASP License Manager" from your computer desktop or double clicking on "lmsetup.exe" in "C:\Program Files\SR Research\Common" folder. Failing to install this will report a "NO HASP Key Found" error in the SR License Manager dialog box.
 - When installing the license manager, set the "Installation Type" to Service (nhsrvice.exe).
 - On the "HASP License Manager" dialog box, check either Yes or No to continue.
4. If this is the server computer, you may install optional Aladin Monitor software by clicking "Start -> All Programs -> SR Research-> Aladdin Monitor" from your computer desktop or double clicking on "Aksmon32.exe" in "C:\Program Files\SR Research\Common" folder.
5. Test license status.
 - Please make sure that the server and client computers are running and visible to each other in the same network group (check this out from "My Network Places -> View Network computers"). Contact your system administrator if the computers cannot see each other in the same network group.

- Please make sure that the network license dongle is plugged to server computer (remove all other HASP dongles) and drivers are already installed.
- Now, click "Start -> All Programs -> SR Research -> License Manager" to check for the licensing status for each of the client computers in the network.

3 Working with Files

The EyeLink Data Viewer can import EyeLink EDF files. The viewer loads each data file into a Data Viewing Session (DVS). A DVS can consist of many EyeLink data files. Only one DVS can be open within the Data Viewer at a time.


Once a DVS has been created and EyeLink data files have been imported into it, the DVS file can be saved and then reloaded at a later date for future analysis. It is *much* quicker to load a DVS than to load the original EDF files of which the DVS consists.

All changes made during an viewing session only affect the DVS data; the original EDF files are never altered.

3.1 Creating a Data Viewing Session (DVS)

From the menu, choose:

File → New


Tip: A new viewing session can also be created by clicking the **Create New Viewing Session** tool  on the standard toolbar, or press **Ctrl N** on Windows or **⌘ N** on Mac OSX.

Tip: To find out the meaning of a tool on the standard toolbar, simply place the mouse over the tool.

3.2 Saving a Viewing Session

From the menus, choose:

File → Save

Tip: The viewing session can also be saved by clicking the **Save Viewing Session** tool  on the standard toolbar, or press **Ctrl S** on Windows or **⌘ S** on Mac OSX.

When a new viewing session is saved, Data Viewer will create a {session name}.evs file as well as a {session}.res folder; both of them are required for a proper re-opening of the viewing session later on.


3.3 Opening an Existing Viewing Session

To open an existing viewing session

- 1) From the menus, choose:

File → Open

- 2) In the Open dialog box, browse to the location of the file and select it.
- 3) Click Open


Tip: A saved viewing session can also be opened by clicking the Open Existing Viewer Session tool  on the standard toolbar, or press Ctrl O on Windows or ⌘ O on Mac OSX.

The viewing session can also be opened by double clicking on the {session name}.evs file. Please make sure that the {session}.res folder is available; otherwise the viewing session cannot be properly opened.

3.4 Saving an Existing Viewing Session to a Different File

To save a viewing session to a file

- 1) From the menus, choose:
File → Save AS
- 2) In the Save As dialog box, browse to the directory in which you want to save the document.
- 3) In the File name text box, type in the name for the document.
- 4) Click Save.


Tip: The viewing session can also be saved by clicking the Save Viewing Session As tool  on the standard toolbar.

3.5 Importing EyeLink Data Files

EyeLink EDF files can be imported into an already opened Viewing session or, if no Viewing session is open, a new one will be created for you when the data files are imported.


To import an .EDF file,

- 1) From the menus, choose:
File → Import Data → EyeLink Data File
- 2) In the Load dialog box, browse to the location of the EDF file that you want to import and select it.
- 3) Click Load.

Tip: EyeLink data can also be imported by clicking the Import EyeLink Data File button  on the standard toolbar,

Multiple EDF files can be imported through “File → Import Data → Multiple EyeLink Data Files”, or Press Ctrl I on Windows or ⌘ I on Mac OSX. In the “Import EyeLink Data File” dialog box, browse to the intended “Location”. A list of EDF files in the

folder and sub-directories will be displayed. Select the intended files and then click “Import”.

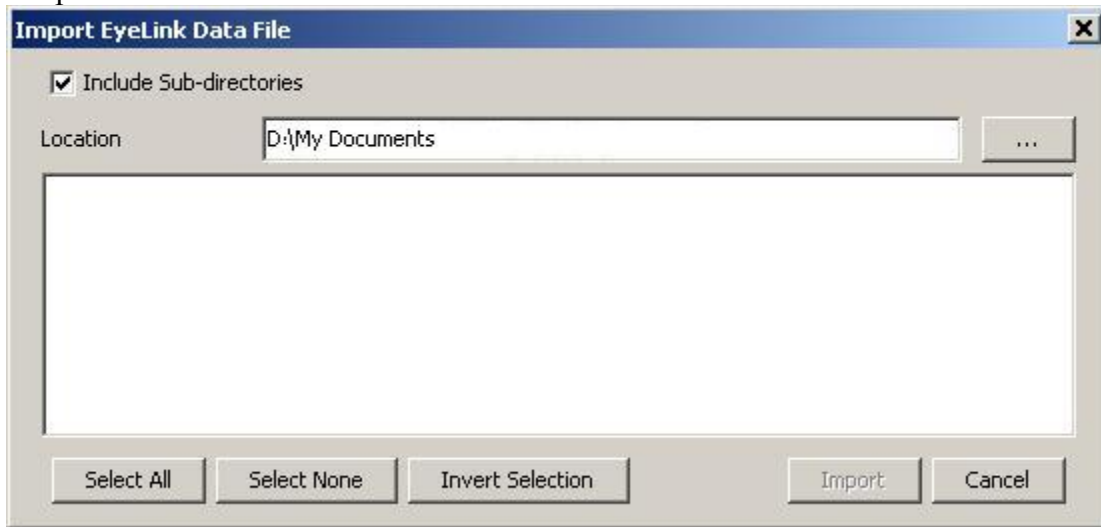

Version 1.7 of Data Viewer introduced a change to optimize the viewer performance by not loading the samples by default. To make sample data available to a viewing session, you will need to first change the data loading preferences and then reload the EDF file.

1. Before loading any EDF files, please close the current viewing session and click "File -> New" to create a new session.
2. Click "Preferences" tab and go to "Data Loading" preferences. Check the "Load Samples" option. If you want this change to be persistent over future viewing sessions, you may consider saving this as default setting. This can be done by selecting the topmost preference tree node, clicking the right mouse buttons and selecting "Save Properties as Defaults".
3. Now start loading in the EDF files.

Note: The time it takes to load the data files will depend on their size and heavily on whether or not you are loading samples from the data file. A 60 minute, binocular, 500Hz recording will take 3 – 4 minutes to initially parse the data file if loading all events and samples. Once you have saved the viewing session, reloading the session takes a fraction of the time because the data is in a format optimized for the Data Viewer.

Note: To load data files using the SceneLink gaze mapping, please make sure to enable the “Enable SceneLink Gaze Mapping” and “Generate SceneLink Frame Messages” in the “Data Loading” preferences settings before you import the data files. Make sure that you have already created the lookup tables for the trials in SceneLink.

4 Data Viewer Windows

The EyeLink Data Viewer uses a desktop framework that contains all windows of the Viewer application. Once a Data Viewing Session (DVS) is opened and EyeLink data is loaded, two windows will appear: the Inspector Window on the left and the Trial View Window on the right. The Inspector Window enables the user to choose data to be viewed, to review statistical summaries, and to configure default application settings. The Trial View window allows a graphic examination of the data of the selected trial. Both windows can be minimized and maximized within the Application desktop. The figure below shows the viewer desktop after loading the sample picture.edf file.

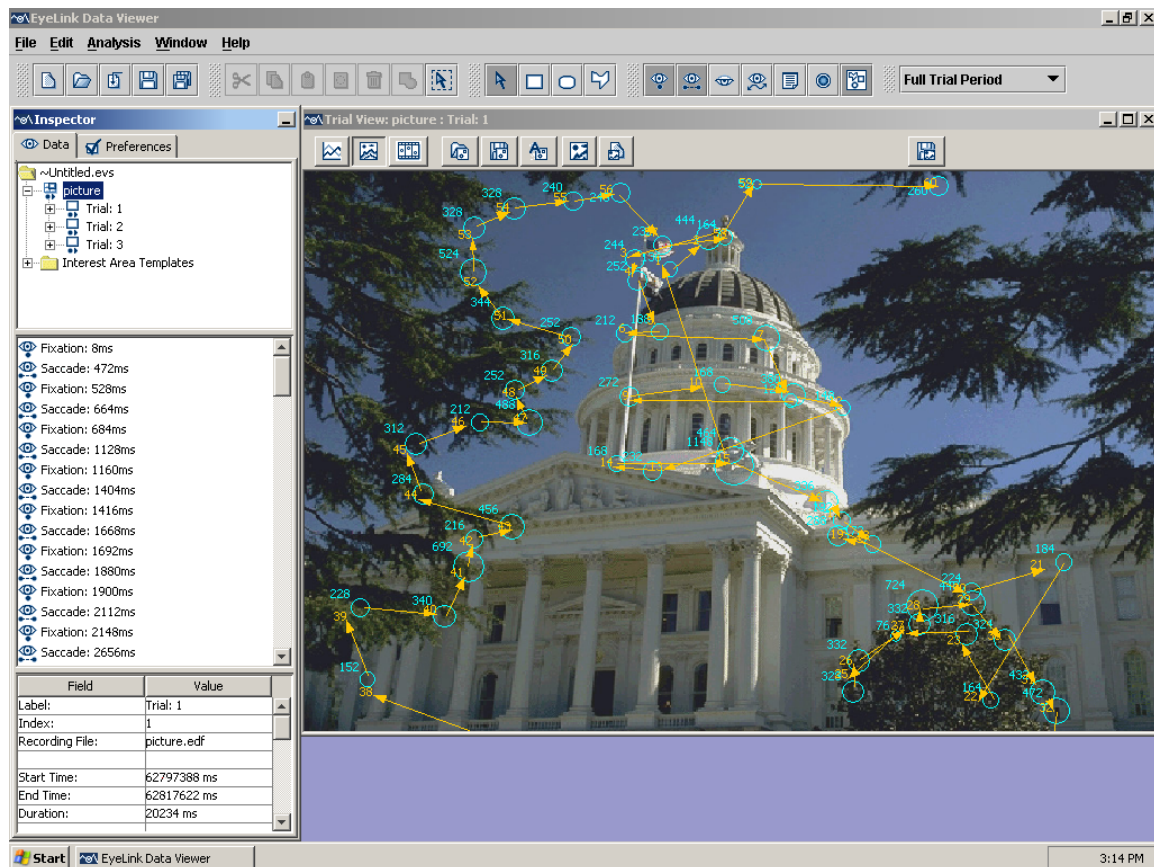


Figure 4-1. Viewer Display after Loading the Sample Picture.edf File

4.1 Inspector Window

The Inspector Window consists of 2 tabs, the Data Tab and the Preferences Tab (see Figure 4-2).

4.1.1 Data Tab

The Data Tab consists of 3 panels (Figure 4-2, left Panel). The top panel provides a tree view of the viewing session. The first level of the tree view contains trial-grouping nodes (). EyeLink Data Viewer lists trials within a viewing session by either data file or by experimental conditions. Each trial-grouping node has a set of tree leaves representing the trials () that are contained within that node.

The middle panel represents a list of components that are within the selected node of the analyses session tree. If a trial is selected in the viewing session tree, a list of visible events is displayed. If an Interest Area Set is selected, a list of interest areas that make up the set are selected.

The bottom panel is a table representing the properties of the selected tree node or selected list component. For example, if a fixation is selected, all properties of the fixation are listed in the properties table. Some properties, like color, are editable, by pressing on the property value with the mouse.

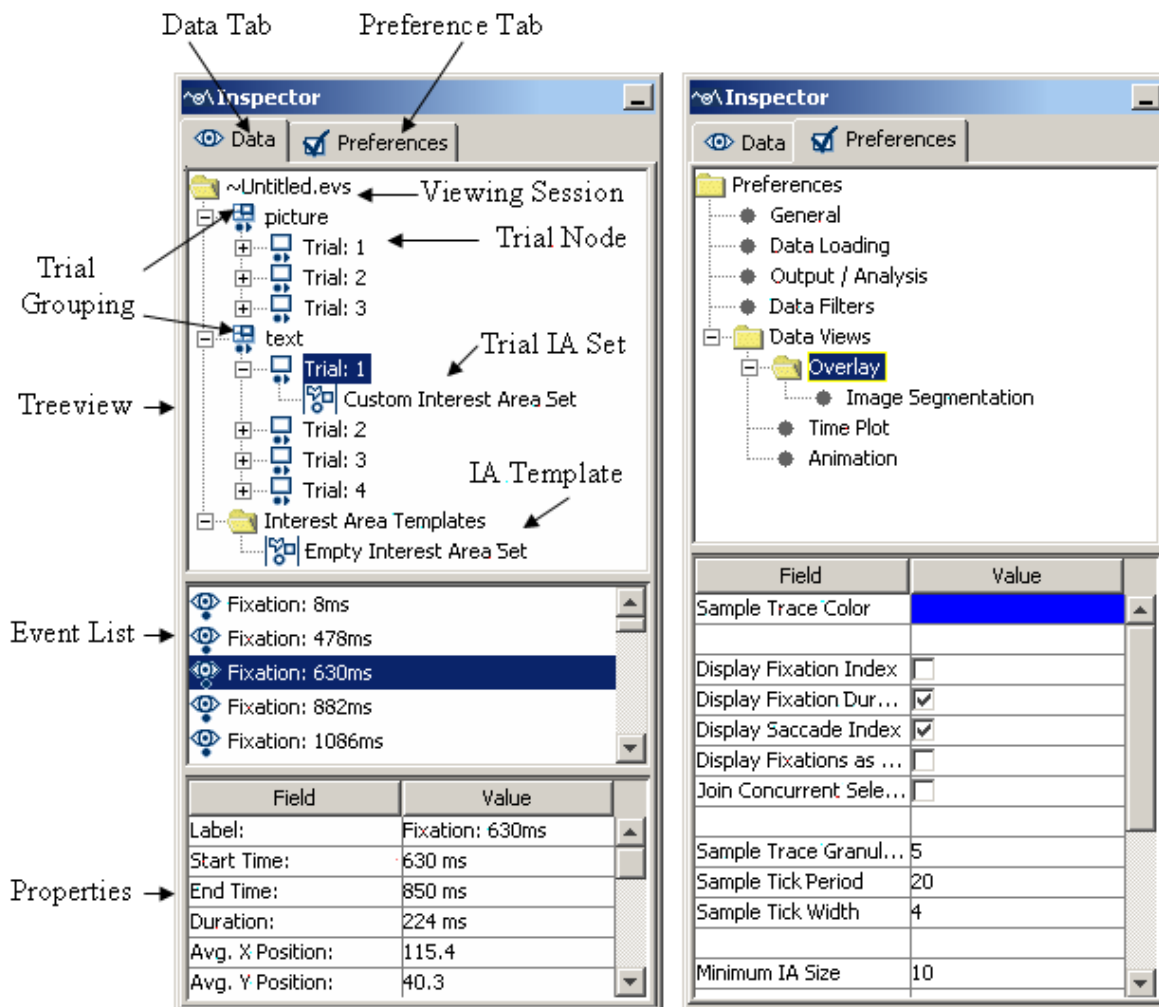



Figure 4-2. Components of the Inspector Window (Left: Data Tab; Right: Preference Tab)

4.1.1.1 Data Tree Hierarchy

When a data file is loaded, the tree view panel lists the possible analysis levels in a hierarchical order. The user can obtain information on the trial groups and individual trials. Within each trial, statistics on all events (fixations, saccades, blinks, button presses, and messages) and interest areas can also be retrieved.

4.1.1.2 Trial Grouping

By clicking on a trial-grouping node () , information on this trial group can be obtained in the bottom panel of the data inspector. Typically, this includes the label of group (values of the grouping variables), total number of trials, and configurable interest area information (see section 6.1.3 on how to perform trial grouping). By default, trials in a viewing session are grouped by the EyeLink data files loaded. The following table lists the group properties if trials are grouped by data file. Those fields that are editable in the viewer are marked with an asterisk.


Field	Contents
Label *	The label of the trial group
Number Trial	The total number of trials in the trial group
File	The file name of the recording (with the path information)
Eyelink Version	Tracker version of the EyeLink system
Sampling Rate	250 or 500 Hz
Eyes	Which eye produced the recording (the value could be Left, Right, or Binocular)
Mode	Pupil only mode, or pupil with corneal reflection
Pupil Data Type	Whether the pupil size is set by area or diameter
Sample Data type	Raw (raw pupil-center position or pupil minus corneal if running in pupil-CR mode as measured by the image-processing system), Gaze (actual gaze position on the display screen, fully compensated for head position), or HREF (the tangent of the rotation angle of the eye relative to the head)
Default Interest Area Set	This shows the default interest area set (chosen from the interest area templates) used for this group of trials. If none, then displays “Empty Interest Area Set”.
Display Width *, Display Height *	Screen resolution (width and height in pixels) of the recording displays. By default, this sets the trial view window resolution to the value specified by display Coordinates messages. If you want to change the screen resolution, make sure that you first

	close the Trial View Window first, change the Display Width and Display Height values, and then open another trial view window.
Report Hidden	If checked, the entire data file will be excluded from the output reports.

4.1.1.2.1 Deleting a group of trials

To remove a group of trials from the analysis, select the grouping node, click the right mouse button and click on the “delete” menu item. If trials contained within a grouping node are deleted from the data viewer session, they will be permanently removed from the tree view (the EDF file has to be reloaded before information on that file can be accessed again).

4.1.1.3 Trial Summary

By clicking on the trial node () of the tree view, information on participants' performance in individual trials can be retrieved. Once a trial is selected for analysis, the middle panel of the Data Tab lists all of the visible events and messages in the trial and the bottom-panel provides a general summary of performance in that trial. The following table lists the trial-level analysis variables. Those fields that are editable in the viewer are marked with an asterisk.

Field	Contents
Label *	Label of the trial
Index	Trial sequence
Recording File	File name of the recording
Recording Start Time	Timestamp when the trial recording started (in milliseconds since EyeLink tracker was activated)
Recording End Time	Timestamp when the trial recording ended (in milliseconds since EyeLink tracker was activated)
Recording Duration	Duration (in milliseconds) of the trial recording
Interest Period Start	Start time of interest period (in millisecond relative to the trial recording start time)
Interest Period End	End time of interest period (in millisecond relative to the trial recording start time)
Reaction Time	Configurable response time determined by the Reaction Time Definition associated with the trial. Use the value of Duration if RT is undefined.
RT Event Index	Index of event that triggered the end of the reaction time; -1

	when NO_MATCH. The index is relative to a specific event type; for example an index of 5 for a Saccadic RT means the 5 th saccade in the trial.
RT Event Type	The way in which the trial reaction time is ended, possible values are TIMEOUT, BUTTON, FIXATION, SACCADE, MESSAGE, or NO_MATCH
RT Definition Label	Label of the reaction time definition (see section 6.3.1 Managing Reaction Time Definitions). “.” if undefined; “CUSTOM” if the RT definition is a custom one (see Section 7.4 Reaction Time Definitions).
Average Fixation duration	Average duration (in milliseconds) of all fixations in the trial
Average Saccadic Amplitude	Average size (in degrees of visual angle) of the saccades in the trial
Report Hidden *	Whether or not to hide the trial from output reports.
Interest Area Set *	Setting of the interest areas (“Custom Interest Area Set” or “Empty Interest Area Set”). If interest areas have been defined, they can be removed by selecting “Empty Interest Area Set”
Number Samples	Total number of samples in the trial recording
Number Fixations	Total number of fixations in the trial recording
Number Saccades	Total number of saccades in the trial recording
Number Blinks	Total number of blinks in the trial recording
Number Button Presses	Total number of button presses in the trial recording
Number Messages	Total number of messages in the trial recording
X Resolution, Y Resolution	Angular resolution at current gaze position (in screen pixels per visual degree)

4.1.1.3.1 Hiding a Trial

To hide a trial from the output reports

- 1) Click the trial to hide in the top panel of the data tab.
- 2) In the bottom panel of the data tab (which contains properties for the trial), find the entry “Report Hidden” and change the current setting.

4.1.1.3.2 Deleting a Trial


To delete a trial from the Data Viewer, simply place the mouse over that trial node, click the right mouse button and select **Delete**. The trial is permanently removed from the data viewing session and all other trial indexes are updated accordingly.

4.1.1.4 Interest Area Templates

An interest area template is a .IAS file that contains a set of interest areas. If the same interest areas are used for several trials, instead of repeating the procedure of creating interest areas for each individual trial, an interest area template can be loaded to supply segmentation information for those trials efficiently. The property panel lists the following summary statistics for an interest area template. Those fields that are editable in the viewer are marked with an asterisk.

Field	Contents
Name *	Label of the custom interest area set
IA Count	Total number of interest areas defined in the trial
IsCustom	Whether the current set of interest areas is user-defined (true) or an interest area template (false)
Image Name Mapping	If the "Auto Select IAS for Trial" property in the "Output/Analysis" preference settings is true, the value in this field will be used to associate the IAS to trials automatically (see section 5.8.8 "Automatic Trial Image-Interest Area Set Association" for details).

4.1.1.5 Custom Interest Area Set/Empty Interest Area Set

For each trial, the tree view also shows whether or not a custom interest area set has been defined for the trial. If there is no interest area set, the node  will be "Empty Interest Area Set". The property panel lists the same elements as the interest area templates, except that the IsCustom field is "true" for the custom interest area set.

4.1.2 Preferences Tab


The preferences tab (Figure 4-2, right panel) provides access to all the elements of the data viewer that can be configured by the user. These include everything from the default colors used for displaying events, to the elements of an EDF file to load. These elements are broken down into the following five categories: General Preferences, Data Loading Preferences, Output/Analysis Preferences, Data Filters Preferences, General Trial View Preferences, Spatial Overlay View Settings, Temporal Graph View Settings, and Animation Playback Settings.








4.2 Trial View Window

The Trial View window displays the data of a selected trial in a Spatial Overlay Mode, a Temporal Graph Mode, or an Animation Playback Mode. Selecting a trial in the Data Inspector tree view shows the trial's data in a Trial View Window. The user can have

several Trial View Windows open at one time by right clicking on a trial in the Inspector window and selecting “New Window” in the popup menu.

4.2.1 Selecting Data Elements to Display

At anytime event visibility can be toggled on/off within the Trial View Window. For example the  button in the Application Desktop toggles the visibility of Fixation Events.

	Toggle Fixation Event Visibility
	Toggle Saccade Event Visibility
	Toggle Blink Event Visibility
	Toggle Message Event Visibility
	Toggle Button Event Visibility
	Toggle Sample Visibility
	Toggle Interest Area Visibility

4.2.2 Cloning a Trial View

From the menu, choose:

Window → Clone Window.

This allows, for example, both the temporal graph and spatial overlay views to be shown at the same time on the desktop. You can tile the windows together. Note that event selection in one view will automatically cause the same events to be selected in the other view.

4.2.3 Tiling/Cascading Windows

To arrange all open trial view windows, from the menu, select:

Window → Tile

or

Window → Cascade

Figure 4-3 shows an example of tiling three open trial view (spatial overlay, temporal graph, and animation modes) windows.

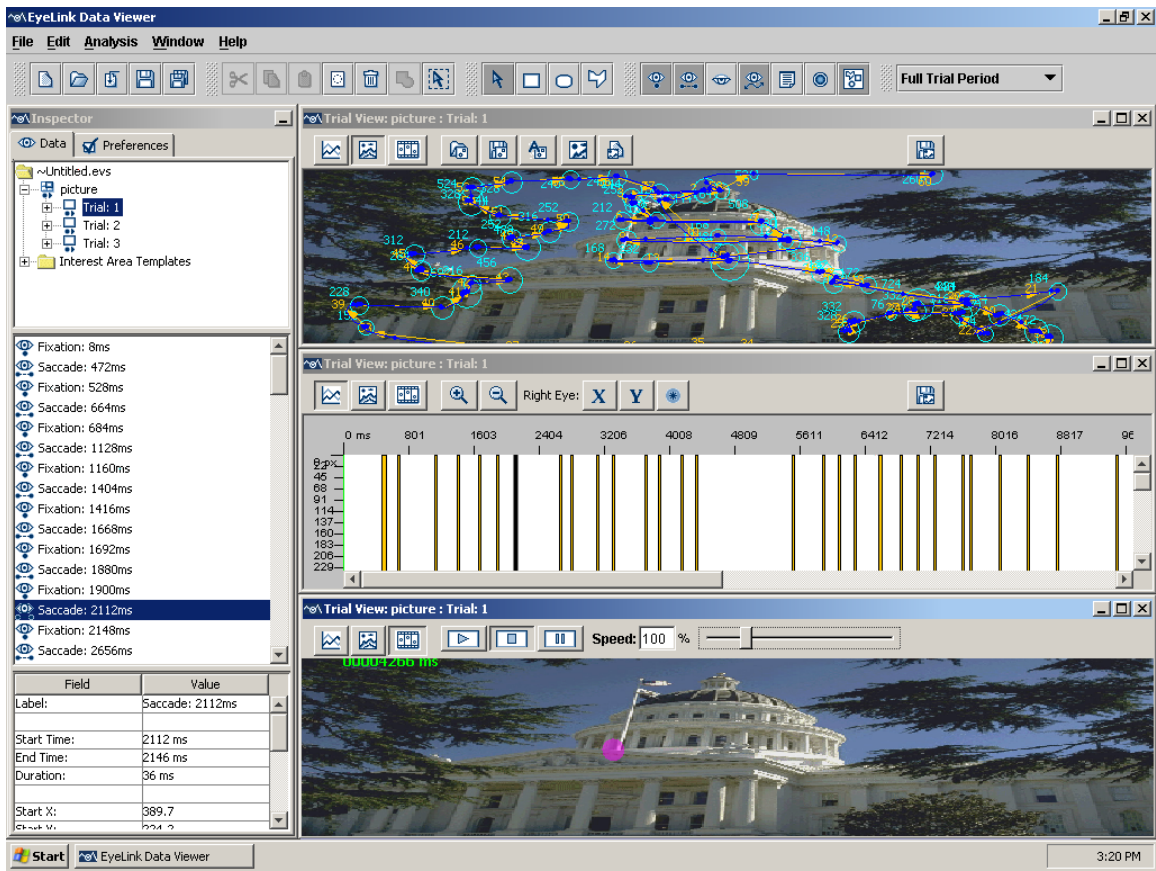



Figure 4-3. Tiling Trial View Windows.

4.3 Spatial Overlay Trial View

The Spatial Overlay view (see Figure 4-4) is selected by pressing the  button in the Trial View window. This view lets you look at events and samples by placing them where they were detected in space. If the image that was presented during the trial is available, it can be placed in the overlay view.

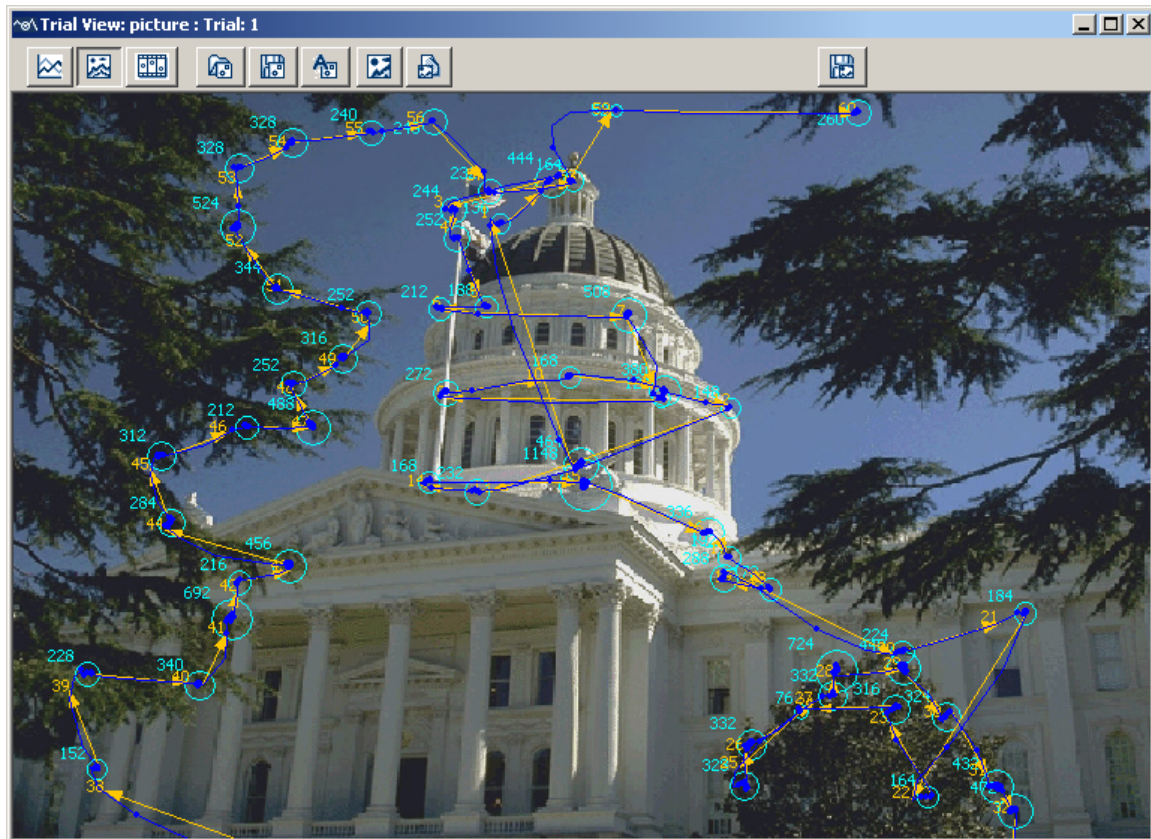



Figure 4-4. Sample Spatial Overlay Trial View

The following buttons are available from the spatial overlay mode:

	View Trial Data in 2D Temporal Graph Mode: Select the temporal graph mode.
	View Trial Play Back Animation: Select the playback animation mode.
	Save Trial View as Image: Creates an image of the spatial overlay view
	Import Area of Interest File: Loads an interest area set file and sets the interest areas for this trial. Any existing interest areas are deleted from the trial. After loading, the Interest Area Set for the trial is set to “Custom”.
	Save the Interest Area Set File: Saves the trials interest area set to a file.
	Auto Segment Image to Interest Areas: Creates interest areas automatically for the trial bitmap.
	Display Image in Full. If enabled, overlay is not scaled to size of Trial Viewer Window and is instead set to actual size (One pixel of overlay corresponds to one pixel of trial display).
	Change the Overlay Image: Select the image to use as the background of the overlay. Any existing background graphics are deleted.

Changes from the previous versions: Note that when the background image for a trial is manually changed with the “Change the Overlay Image” button, please also check for the “Fit Overlay Image Change to Display” setting in the Data Views->Overlay preference. If that setting is true, the image is scaled to fit the display dimensions for the trial. If false, the image size is not altered and the image is centered on the trial overlay view.

4.4 Temporal Graph Trial View

The Temporal Graph View (see Figure 4-5) allows the user to view trial data as a trace plot, where the X axis represents time and the Y axis represents the X / Y gaze location, pupil size (divided by a factor of 10), acceleration, and velocity data. To view trial data in this mode, select the  button in the top left corner of the Trial View Window.

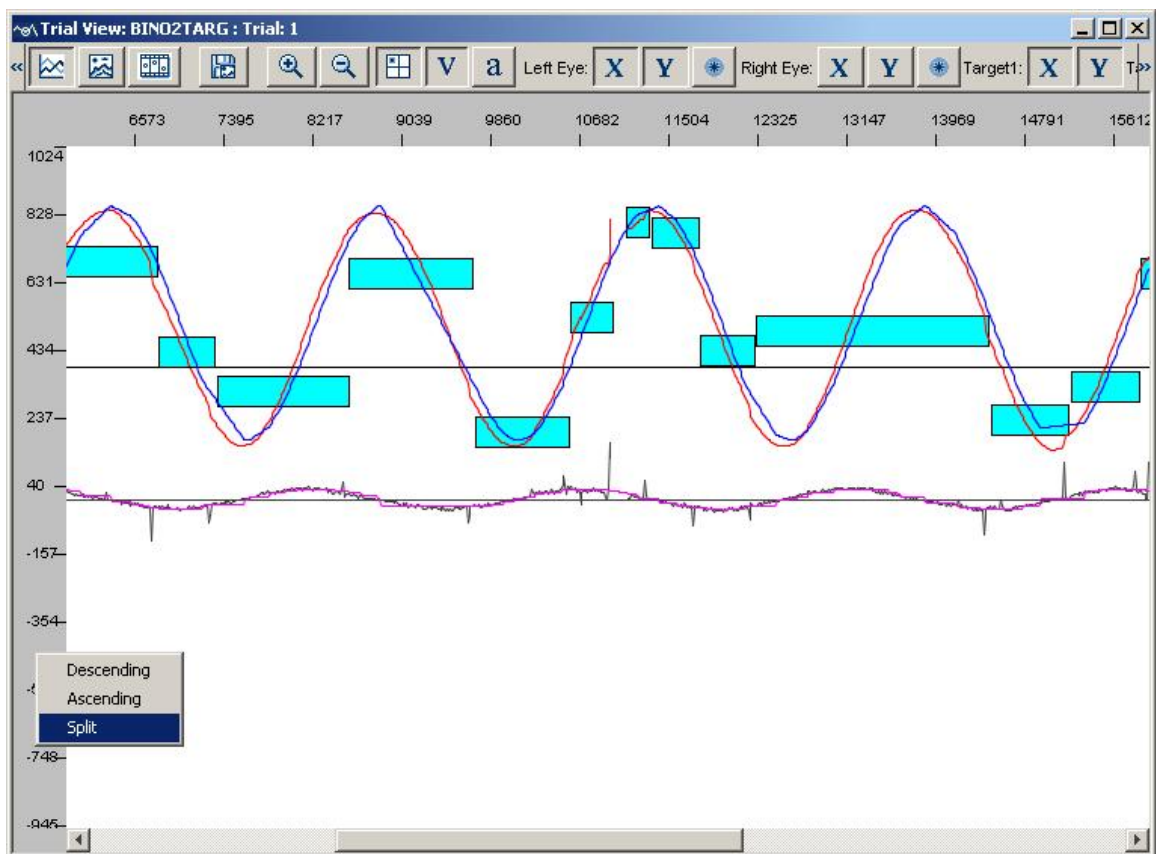














Figure 4-5. Sample Temporal Graph Trial View

4.4.1 Viewing Sample and Target Data

Time plot view will show trace plots only if the EDF file contains samples and if sample are activated for viewing (i.e., the “Toggle Eye Sample Visibility” button  on the Data

Viewer tool bar is selected). The position, velocity, and acceleration sample traces can be selected for viewing from buttons within the Trial Viewer Window.

	View Trial Spatially with Image Overlay: Select the spatial overlay mode
	View Trial Play Back Animation: Select the playback animation mode
	Save Trial View as Image: Creates an image of the temporal graph view
	Zoom In Graph View: Zoom in to view details of a selected portion of data
	Zoom Out Graph: Zoom out the graph by a scale of 2
	Position Trace: Show the gaze or target position.
	Velocity Trace: Show the gaze or target velocity data
	Acceleration Trace: Shows the gaze or target acceleration data
	X Sample Trace: Toggle X trace visibility
	Y Sample Trace: Toggle Y trace Visibility
	Pupil Sample Trace: Toggle Pupil Trace Visibility

Changes from the previous versions: to show the eye or target traces in this version, the user needs to select the type of data (position, velocity, or acceleration) and data source (x, y, pupil of left eye and/or right eye, x, y of target1 and/or target2) to be shown. To show the position, velocity, and acceleration traces of target(s) in this view, messages on the current position of the target(s) must be recorded in the EDF file. These messages should contain a target position keyword (“!V TARGET_POS” by default and can be configured in the “target position string” of Data Loading preference settings), followed by parameters of the individual targets.


The scale in which the vertical data is displayed can be changed by clicking the right mouse button when the mouse cursor is placed on the Y-axis and choose one of the following options: ascending (large values on the Y-axis appear on the top of the screen), descending (large values on the Y-axis appear at the bottom of the screen), or split (0 appear in the middle of the screen, positive values on the top and negative values at the bottom).

The X, Y, and pupil traces can be highlighted. Press the 'x' key to highlight the X trace, press the 'y' key to highlight the Y trace, and press the 'p' key to highlight the Pupil trace. If pressing one of these keys does not highlight the intended sample traces, please make sure that the trial view window is currently in focus (double clicking anywhere the sample traces are plotted to set the keyboard focus to the trial view window).


Various aspects of the time plot view can be configured in the preference settings (Preferences -> Data Views -> Time Plot). For example, the user may change the colors of the sample traces, the scaling factor for plotting the acceleration, velocity, and pupil size data, total number of initially visible samples in the temporal graph view window, etc.

4.4.2 Zooming Operations


The temporal graph can be zoomed in and out. To zoom into an area of the temporal graph:

- 1) Hold the Shift key down or press the  button.
- 2) Position the mouse on the temporal graph at the location of the upper left hand corner of the area you wish to zoom into.
- 3) Press the mouse button and drag the mouse to the lower right hand corner of the area you wish to zoom into.
- 4) Release the mouse button
- 5) Release the Shift key.








The graph will be magnified so that the area you selected fills the Trial View Window.

To zoom out the graph, press the  button. The temporal graph will be zoomed out by a factor of 2.

4.5 Animation View

The Animation View (see Figure 4-6) is selected by pressing the  button in the Trial View window. This view plays back the gaze position (superimposed on the background image that the participant was looking at while the data was being recorded), with a concurrent time code displaying the time from the start of trial recording.

Several controls are available for the animation view. The user can start, stop, or pause the playback by clicking the appropriate button. The speed of playback can also be adjusted (the user modifies the speed value and then press the “Enter” key to register the change). A trackbar control displays the progress of the playback and lets the user select a position to start playback.

	View Trial Spatially with Image Overlay: Select the spatial overlay mode
	View Trial Data in 2D Temporal Graph Mode: Select the temporal graph mode
	Playback Trial: Starts playing back the trial data
	Stop Trial Playback: Stops the playback and resets the time code and gaze cursors.
	Pause Trial Playback: Pauses the playback
	Forward Step: Each click on this button advances the playback by one video frame.
	Save Trial to Movie File: Save the playback of the trial into a movie file

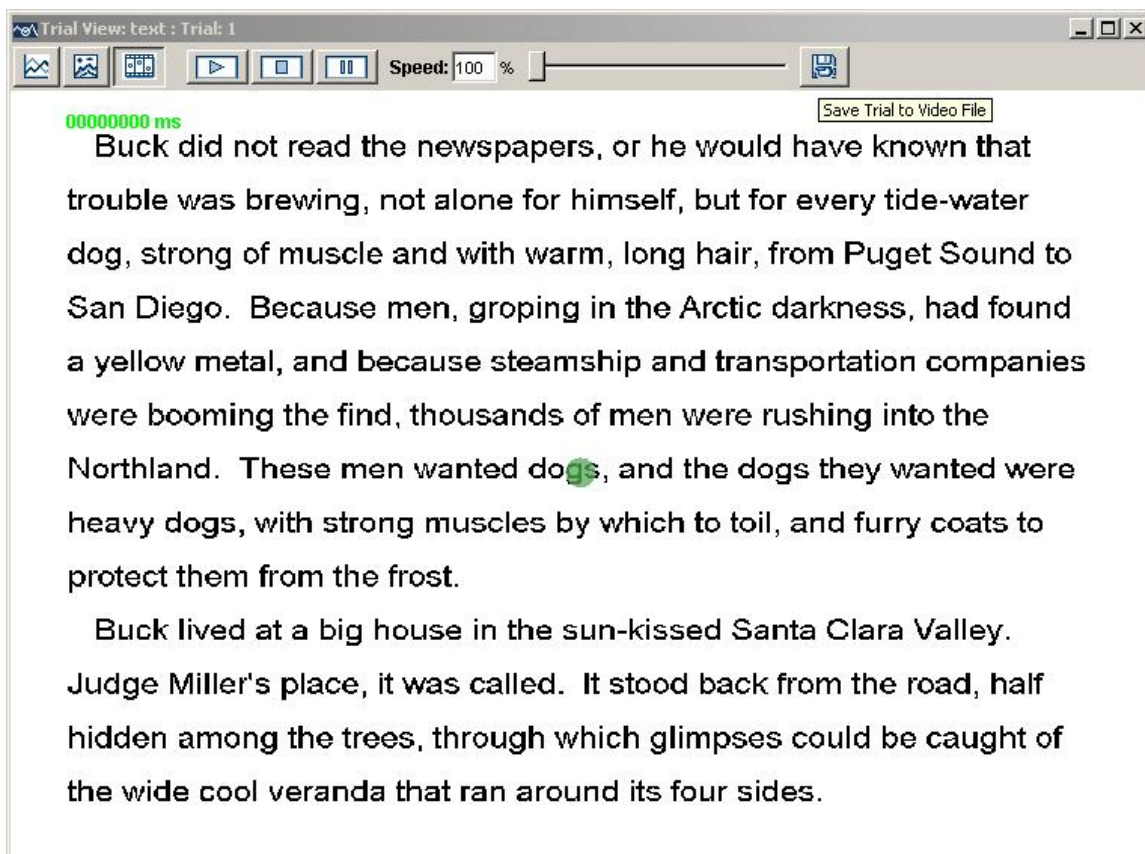



Figure 4-6. Sample Playback Animation View.

To save the playback of trial data into a movie file, click on the "save To Movie" button () on the trial view window toolbar. In the following "Movie Save Options" dialog box, enter the movie file name, file type (.AVI or .MOV), frame rate, and compression method. Click "Start" to begin moving saving. Depending on the size of recording file, this may take a long time to finish. At any point, user can press "ESC" key to stop the saving process.

Playing back video clips in the animation view is also supported if the EDF file is created from an SR Research Experiment Builder recording. The video clip must be encoded in XVID format and a computer with P4 CPU is required for playback in Data Viewer. The current frame number is displayed at a position relative to the top-left corner of the video. If you want to hide this frame number indicator somehow, please go to "Preference Settings -> Animation" and uncheck the "Display Frame Number" option. Please be advised that this feature is presently not available on Mac OSX.

If you want to save the video clip with the overlaid eye position, please try to make sure the display solution on your computer desktop (right click on the desktop to start the "Display Properties") matches or is close to that of the EDF file recording resolution (the "Display Width" and "Display Height" in the recording session node) for best video

saving quality. In the video saving dialog box, set the "size" to this recording resolution. The default NONE compression will save a large output file. You may try some codec that does the compression (e.g., the "CINEPAK CODEC", which does a good compression albeit slow). Please also keep the default Trial View Window size as any change in the aspect ratio of this window may introduce distortion to the saved overlay video. When the video saving is in progress, please do not disrupt it (e.g., by resizing or moving the Data Viewer window or have it occluded) as whatever is visible on the display frame will be saved in the capture as well.

5 Working with Events, Samples, and Interest Areas

The EyeLink Data File (EDF file) contains two streams of data: eye-position samples (up to 2000 per second produced from the EyeLink tracker, depending on the system model) and events (eye-movement events such as saccades and fixations, subject responses, and synchronizing events from the experimental application). Both streams are time-synchronized for easy analysis. The file is organized into trials of data, one for each of the start and stop of tracker recording. Each trial may have samples, events, or both.

Samples are time-stamped in milliseconds and contain monocular or binocular eye-position data - for now Data Viewer only report the eye position in screen coordinate (GAZE) while other data formats (PUPIL or HREF) are ignored. Pupil sizes as area or diameter are also recordable. Samples may also contain eye-movement resolution (used to compute true velocity or saccadic amplitudes).

Eye-movement events record eye position changes identified by the EyeLink tracker's on-line parser, such as fixations, blinks, and saccades. Both the onset and end of these events are marked, allowing samples to be assigned to eye-movement periods without complex algorithms. Important data for analysis such as average position for fixations and peak velocity for saccades is also recorded in the events. Other events record subject responses (such as button presses) and synchronization and data messages from applications. These can be used to record the time of a change in the display, or an experimental condition.

Events are generated by the EyeLink eye tracker (EyeLink I, EyeLink II, and EyeLink CL) in real-time based on an internal heuristic saccade detector (using a velocity and acceleration-based saccade detection method). A blink is defined as a period of saccade-detector activity with the pupil data missing for three or more samples in a sequence. A fixation event is defined as any period that is not a blink or saccade. To detect a saccade, for each data sample, the parser computes instantaneous velocity and acceleration and compares these to the velocity and acceleration thresholds. If either is above threshold, a saccade signal is generated. The parser will check that the saccade signal is on or off for a critical time before deciding that a saccade has begun or ended. A detailed discussion on event parsing can be found in section "4. Data Files" of the EyeLink II or EyeLink 1000 User Manual.


In addition to providing a file-based and a trial-based summary of performance, the EyeLink Data Viewer allows a detailed examination of event data such as fixations, saccades, blinks, buttons, messages, as well as raw sample data.

5.1 Common Operations

The EyeLink Data Viewer allows the user to carry out various manipulations over the events, such as selecting a portion of data for further scrutiny, excluding or deleting a subset of events from analysis, performing drift correction on fixations, merging neighboring fixations and saccades, etc. Selecting, hiding, and deleting operations are common to all event types (fixations, saccades, buttons, blink events, and messages) whereas merging and drift correction is available to fixation and saccade events only. The common operations are discussed in this section whereas the unique operations pertaining to certain types of events are documented in the relevant event section.

5.1.1 Selecting Events


To select an event (fixation, saccade, blink, button, or message), place the mouse cursor over the event either in the spatial overlay/temporal graph view or in the event list of the inspector window and click the left mouse button.

To select more than one event, hold down the CTRL key on Windows or the command key on Mac OSX and select the target event, until all desired events are selected. If all events in a trial are to be selected, simply click the **Select All Visible Item** button  on the toolbar (or click the right mouse button and select “Select All”).


To select a group of events within a square region, place the mouse cursor at the top-left corner of the region. Click and hold down the left mouse button, drag the mouse cursor to the bottom-right corner of the square region, and then release the left mouse button.

Tip: Holding the CTRL or command key while selecting an event leaves previously selected events selected. If the CTRL or command key is not pressed, previously selected events are unselected.

5.1.2 Deleting Events

Select the events and click the **Delete Selected Items** button  on the toolbar (or click the right mouse button and select “delete”). If an event is deleted, it will be permanently removed from the temporal graph view, spatial overlay view, as well as the output reports, and cannot be “undeleted”. In addition, deleting a fixation or saccade event will have an impact on other temporally contiguous events. For example, deleting a fixation will merge the saccades on either side of the fixation into one saccade. Similarly, deleting one saccade will merge the fixations on either side of the saccade into one big fixation.

5.2 Fixations

To examine the fixation information in the data file, first make sure that the fixation events are visible. This is done by ensuring that the Toggle Fixation Event Visibility button  in the application desktop toolbar is selected.

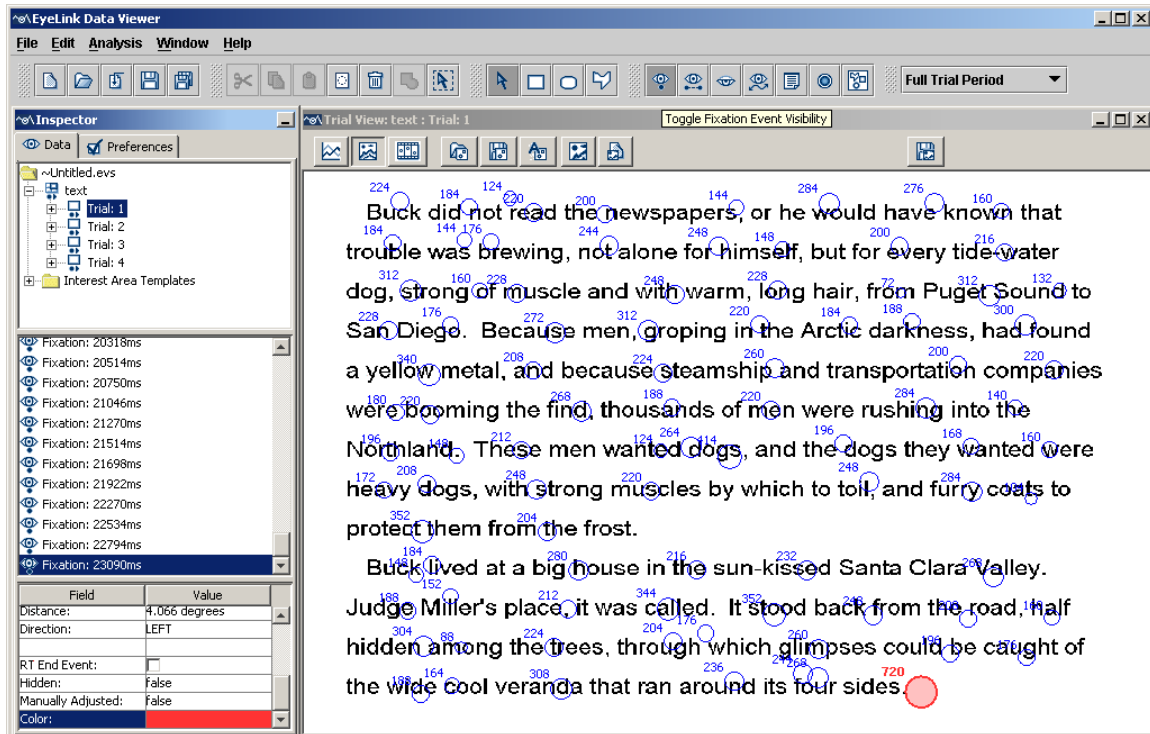


Figure 5-1. Sample Spatial Overlay View of Fixation Events

The trial-view window (Figure 5-1, right panel) shows either a spatial overlay view or a temporal graph view of the fixation events. In the former view, fixations are represented as a hollow circle in the corresponding x, y coordinates on the trial bitmap, with fixation duration displayed near the top-left corner of the circle. In the temporal graph view, fixations are represented as a filled block. All fixation events in a selected trial are listed sequentially in the middle panel of the inspector window. Properties of a selected fixation are shown in the bottom panel of the inspector window. Each field of the fixation event property is described in the following table. Those fields that are editable in the viewer are marked with an asterisk.

Field	Contents
Label *	Label of the fixation
Eye	The eye to which the current fixation event is associated.
Start Time	Trial time ¹ when the fixation starts. Please note that if an interest period filter is applied, the duration, start time, and end time of the fixations that overlap with the start or end of the interest period will be adjusted accordingly.

End Time	Trial time ¹ when the fixation ends.
Duration	Duration (in milliseconds) of the fixation. The fixation duration may be adjusted if the fixation spans the start or end of an interest period and the "Trim Spanned Fixation Time" option is enabled in the Data Filter preferences.
Avg. X Position *, Avg. Y Position *	Average camera X, Y of pupil center
Angle	Angle between the horizontal plane and the line connecting the current fixation and the previous/next fixation
Distance	Distance (in degrees of visual angle) between the current fixation and the previous/next fixation
Direction	Direction ("LEFT", "RIGHT", "UP", or "DOWN") ³ relative to the current fixation in which the previous/next fixation is located
RT End Event *	Whether the current fixation is the reaction time end event (see section 6.3.1 for reaction time definition)
Hidden	Whether the current fixation is hidden from viewing and output analysis. This feature was disabled since version 1.6.
Manually Adjusted	Whether the fixation has been shifted manually
Color *	To achieve better visibility, the color of the currently selected fixation can be changed by clicking on the color toolbox

Note ¹ : System time refers to the timestamp (in milliseconds) since the EyeLink tracker was activated whereas the trial time refers to the time since the trial started. The default trial start time is the start of the trial recording (i.e., "START" message). If reaction-time definitions are applied, the trial-start time is reset at the start event of the RT definition (e.g., "SYNCTIME" message). A negative trial time means that the event or message occurred before the trial start time.

Note ² : The calculation of the angle between the previous/next fixation and the current fixation is directional, with the arrow pointing towards the previous/next fixation (see Figure 5-2 for an illustration of angle calculation).

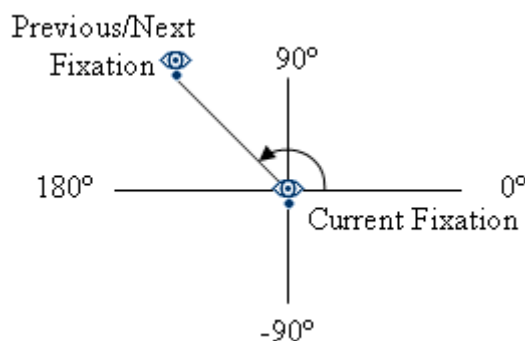



Figure 5-2. Diagram for Angle Calculation

Note³: The direction of the previous/next fixation relative to the current fixation is determined as “LEFT” if the angle is greater than 135° or less than -135°, “UP” if the angle is between 45° and 135°, “RIGHT” if the angle is between -45° and 45°, and “DOWN” if the angle is between -135° and -45°.

5.2.1 Selecting Fixations

In addition to the common method of event selection by mouse, a group of temporally contiguous fixations may also be selected by using a combination of the Shift key and the up or down cursor keys. Select the first fixation by mouse in the event list of the inspector window and hold down the Shift key. Press the down cursor key to select the next fixation, until all target fixations are selected. Use the up cursor key to unselect the recently selected fixations or go to fixations before the first selection.

5.2.2 Merging Fixations

In some cases, two or more temporally contiguous fixations can be merged and represented as a “larger” fixation. To merge fixations, select the fixations and click the Merge Selected Items button  on the toolbar (or click the right mouse button and select “merge”).

Note: Only two or more temporally contiguous fixations can be merged. If fixations are selected that are not temporally contiguous, only the neighboring fixations will be merged.

5.2.3 Drift Correcting Fixations

Sometimes, a drift in the gaze position may occur due to poor system setup, headband slippage, or excessive head/body movement. Normally, the eye tracker operator would notice this problem and perform a recalibration or drift correction. If however, the issue was not addressed during recording, it may be possible to manually adjust the fixation events so that some degree of data analysis can still be performed.

5.2.3.1 Drift-correcting a Group of Selected Fixations

To perform drift correction:

- 1) Select the fixations to be drift corrected
- 2) Hold the ALT key down
- 3) Click anywhere on the trial view window so that it receives the input focus
- 4) Move the selected fixations with the cursor keys to the desired locations

Tip: In case several fixations on one single line (such as in reading research) are to be drift-corrected, another method is to select all of the desired fixations and press CTRL+D on Windows (or Command D on Mac OSX) to align them to the same Y position (see

Figure 5-3). This operation will fail if one or more fixations deviate more than 30 pixels in the Y dimension from the mean of the group of fixations to be drift-corrected in a batch (for example when fixations from different lines are selected). The tolerable deviation threshold and the X/Y dimension of alignment is set to 30 as a default and can be modified in the general trial view preference settings.

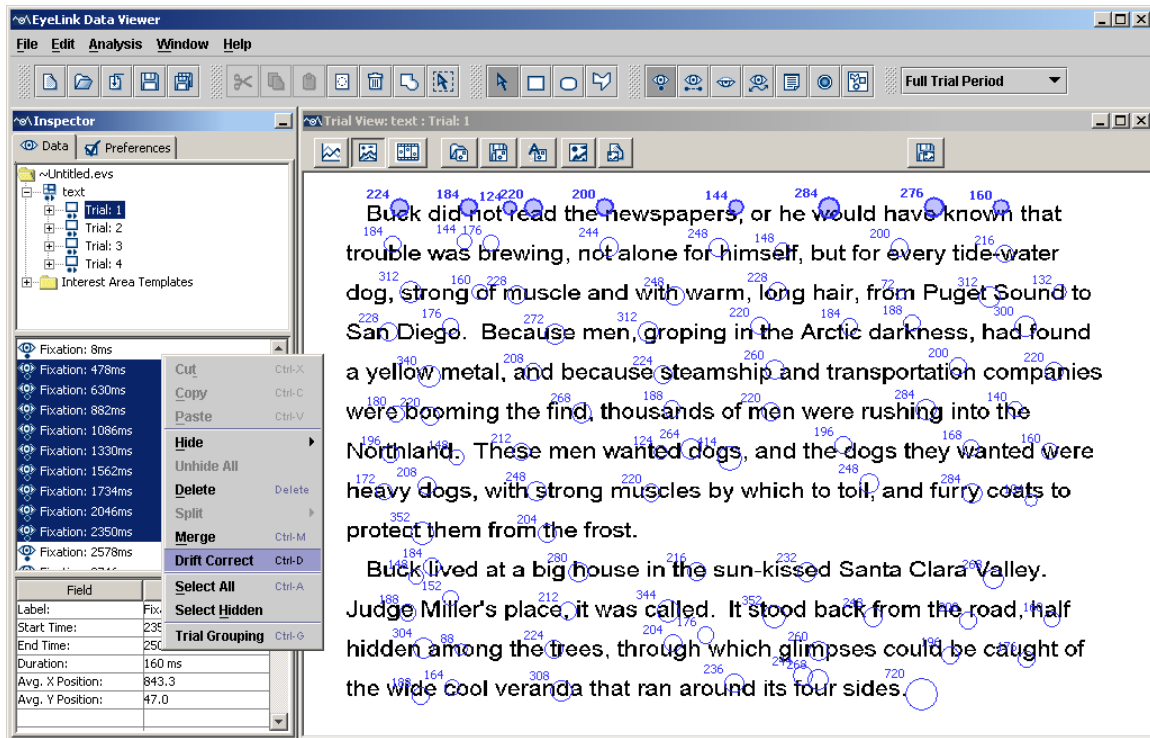



Figure 5-3. Drift Correcting Fixations on One Line of Text.

5.2.3.2 Precise Drift Correction of a Single Fixation

To carry out more precise drift correction of one fixation,

- 1) Select the fixation to be drift corrected
- 2) In the property panel of the Inspector window, select the entry for “Avg. X Position” and type in the desired X position
- 3) Select the entry for “Avg. Y Position” and enter the desired Y position

Note: Drift correcting one fixation or a group of fixations changes the properties of some relevant events, such as the start/end point of a preceding or following saccade and its amplitude.

Tip: Before moving the fixations, it is desirable to turn off the interest area set by pressing Toggle Interest Area Visibility button  in the application desktop toolbar. Otherwise, the interest area(s) within which the fixations appear will also be moved along.

5.2.3.3 Cleaning Fixations

For some studies, researcher may want to clean up fixations based on certain rules. Data Viewer provides two options for fixation cleaning.

1) *Fixation Filter preferences*

The user can filter out small fixations with "Merge Nearby Fixations", "Fixation Merging Amplitude Threshold", and "Fixation Duration Threshold" settings from the Data Filter preference. If the "Merge Nearby Fixations" option is checked, a fixation with a duration shorter than the value set in the "Fixation Duration Threshold" will be merged with a neighboring fixation, if the latter is within the distance set by the "Fixation Merging Amplitude Threshold".

2) *Four-stage Fixation Cleaning (Important: This option was not intended for all users!)*

For some reading studies, the users may want to apply a more refined fixation cleaning algorithm. This can be done by select the intended Trial node or Recording session node, click on the right mouse button, and select the "Clean" option.

	Stage 1	Stage 2	Stage 3
Duration Threshold	80	40	140
Distance Threshold	0.5	1.25	

Stage 4 Durations

Minimum	140
Maximum	800

☐ Delete Fixations Outside Interest Areas

Start

Figure 5-4. Options for Fixation Cleaning.

- i) STAGE 1 For each interest area, Data Viewer checks for each fixation whether its time duration is shorter or equal to the duration threshold. If not, the current fixation will not be cleaned; otherwise, Data Viewer further checks the duration and distance of the fixation immediately before and after the current fixation. The fixation will be merged to the neighboring fixation if its duration is longer than the threshold value and its distance along the x-axis (in degrees) from the current fixation is shorter or equal to the distance threshold. If both the previous and next

- fixations meet the above criterion, the current fixation will be merged to the longer one of the two.
- ii) STAGE 2 This stage is similar to STAGE 1 except for a different fixation duration and a distance value. For this stage to be effective, a shorter duration and a longer distance threshold should be used compared to Stage 1.
 - iii) STAGE 3. Data Viewer goes over the entire data and checks for interest areas which include at least three fixations smaller than the threshold duration and no fixations larger than the threshold duration. In such cases the shorter fixations are merged into the larger ones.
 - iv) STAGE 4 Data Viewer goes over the entire data and deletes every fixation which is shorter than Minimum duration or larger than the maximum duration threshold.
 - v) If "Delete Fixations Outside Interest Area" box is checked, this will remove those fixations falling out side of any interest area. Important: If no interest area is defined, all fixations will be removed!

5.3 Saccades

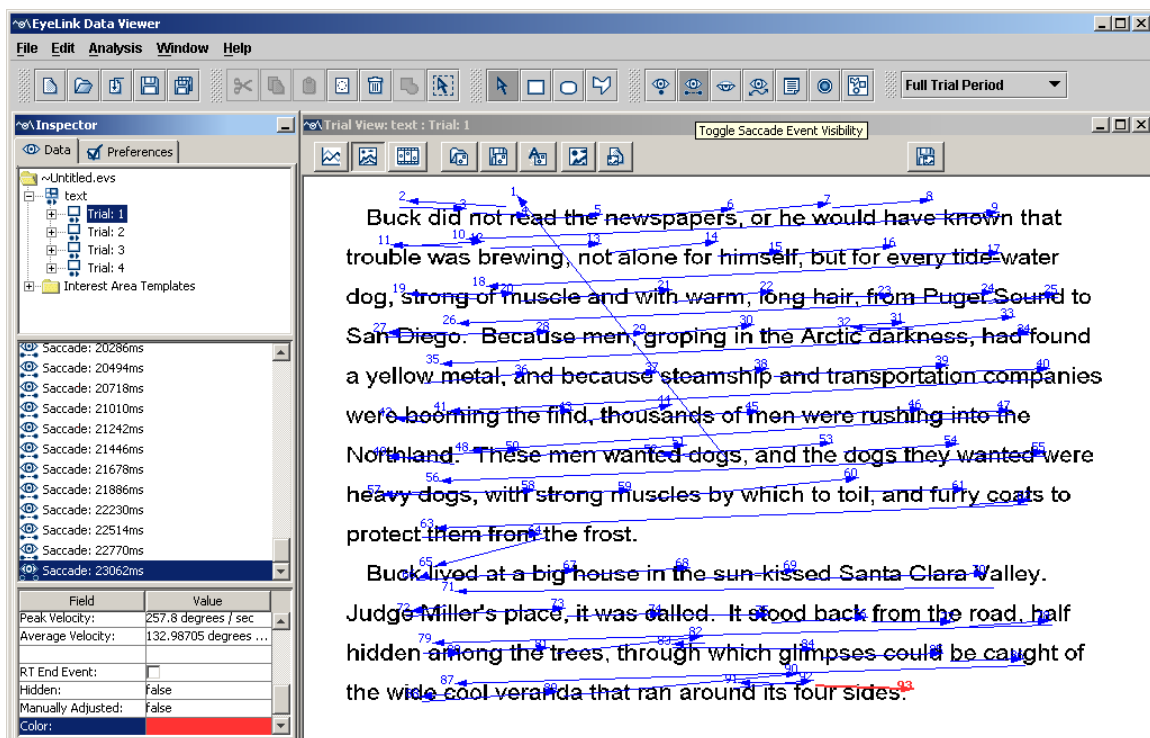




Figure 5-5. Sample Spatial Overlay View of Saccade Events

Saccade events (see Figure 5-4) can be toggled on/off in both the spatial overlay and temporal graph views by clicking the Toggle Saccade Event Visibility button . A saccade is represented as an arrow in the spatial overlay view (pointing towards the endpoint) and as a colored solid line in the temporal graph view. The following table

lists the properties of a saccade event. Those fields that are editable in the viewer are marked with an asterisk.


Field	Contents
Label *	Label of the saccade
Eye	The eye to which the current saccade event is associated.
Start Time	Trial time when the saccade starts
End Time	Trial time when the saccade ends
Duration	Duration (in milliseconds) of the saccade
Start X *, Start Y *	X, Y coordinates of the start position
End X *, End Y *	X, Y coordinates of the end position
Amplitude	Size (in degrees of visual angle) of the saccade
Peak Velocity, Average Velocity	The average and peak values of gaze velocity in visual degrees per second
RT End Event *	Whether the current saccade is the reaction time end event (see section 6.3.1 for reaction time definition)
Hidden	Whether the current saccade is hidden from viewing and output analysis. This feature was disabled since version 1.6.
Manually Adjusted	Whether the saccade has been adjusted manually
Color*	To achieve better visibility, the color of the currently selected saccade can be changed by clicking on the color toolbox.

5.3.1 Merging Saccades

Sometimes, a large saccade is followed by a small corrective saccade or vice versa. Thus, two or more temporally contiguous saccades could be merged. This can be done by selecting the saccades and clicking the **Merge Selected Items** button  on the toolbar (or clicking the right mouse button and selecting “merge”). Once the saccades are merged, information on the ordinal sequence for the current and subsequent saccades is updated and the fixation between the two merged saccades is deleted.

Note: Only two or more temporally contiguous saccades can be merged. If saccades are selected that are not temporally contiguous, only the neighboring saccades will be merged.

5.4 Blink Events

Blink events can be toggled on/off by clicking the **Toggle Blink Event Visibility** button . Blink is represented as a colored solid line in both the spatial overlay view and the temporal graph view. Each field of the blink event properties is listed in the following table. Those fields that are editable in the viewer are marked with an asterisk.

Field	Contents
Label *	Label of the blink
Eye	The eye to which the current blink event is associated.
Start Time	Trial time when the blink starts
End Time	Trial time when the blink ends
Duration	Duration of the blink
Hidden	Whether the current blink is hidden from viewing. This feature was disabled since version 1.6.
Color *	To achieve better visibility, the color of the currently selected blink can be changed by clicking on the color toolbox.

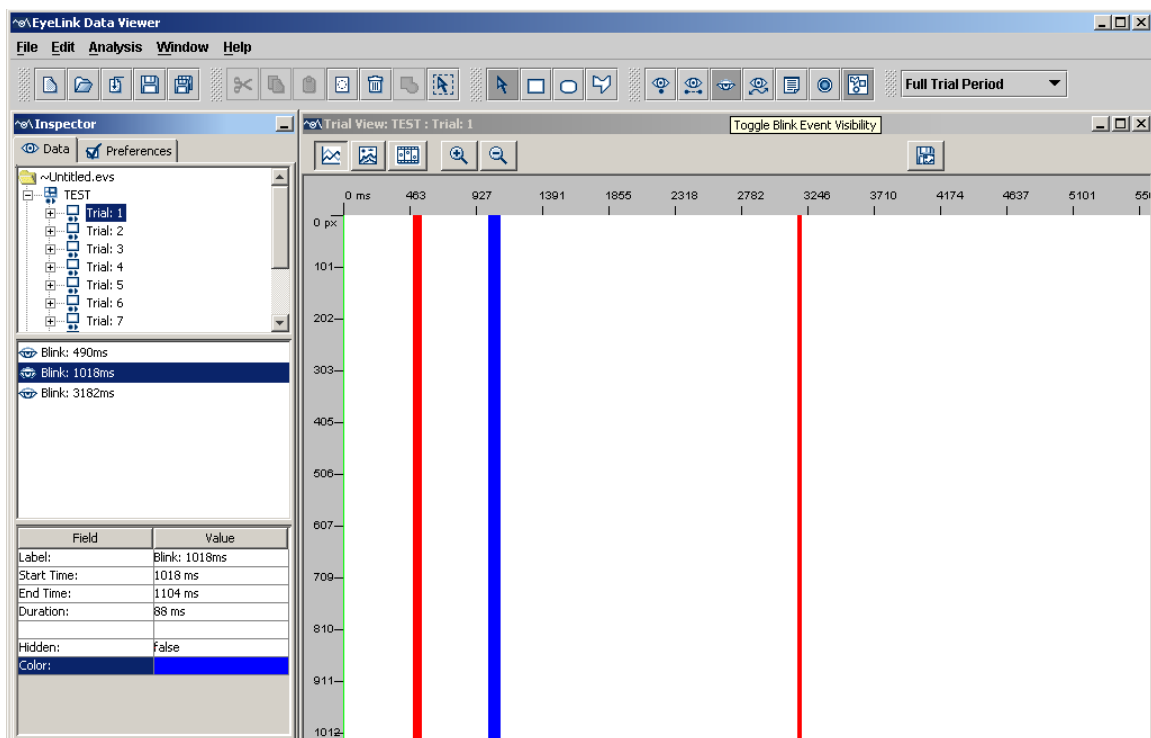



Figure 5-6. Sample Temporal Graph View of Blink Events

5.5 Messages

Messages, including message for built in EyeLink commands, can be placed in the EDF file to mark the start and end of a trial or important events such as display changes. Message events can be toggled on/off by clicking the **Toggle Message Event Visibility** button  in the application desktop toolbar. A message is represented as a colored triangle in the temporal graph view and a colored square in the spatial overlay view, centering on the current gaze position. Each field of the message event property is listed

in the following table. Those fields that are editable in the viewer are marked with an asterisk.

Field	Contents
Label *	Label of the message
Time	Trial time when the message was presented. Note that if the message is presented before the trial starts, a negative time value will be displayed.
Command	If the message is a command message, this lists the actual command used. The commands include simple drawing, built in audio messages, interest area commands, etc. The display commands and audio commands are not shown in the message list unless the “Include Display Command Messages” and “Include Display Command Messages” settings in the Data Filter preferences have been enabled.
Text	Message text
X, Y	Concurrent gaze coordinates when the message is recorded (-100 if no gaze information is available for the time of the message event).
RT End Event *	Whether the current message is the reaction time end event (see reaction time definitions)
Hidden	Whether the current message is hidden from viewing. This feature was disabled since version 1.6.
Color *	To achieve better visibility, the color of the currently selected message can be changed by clicking on the color toolbox.

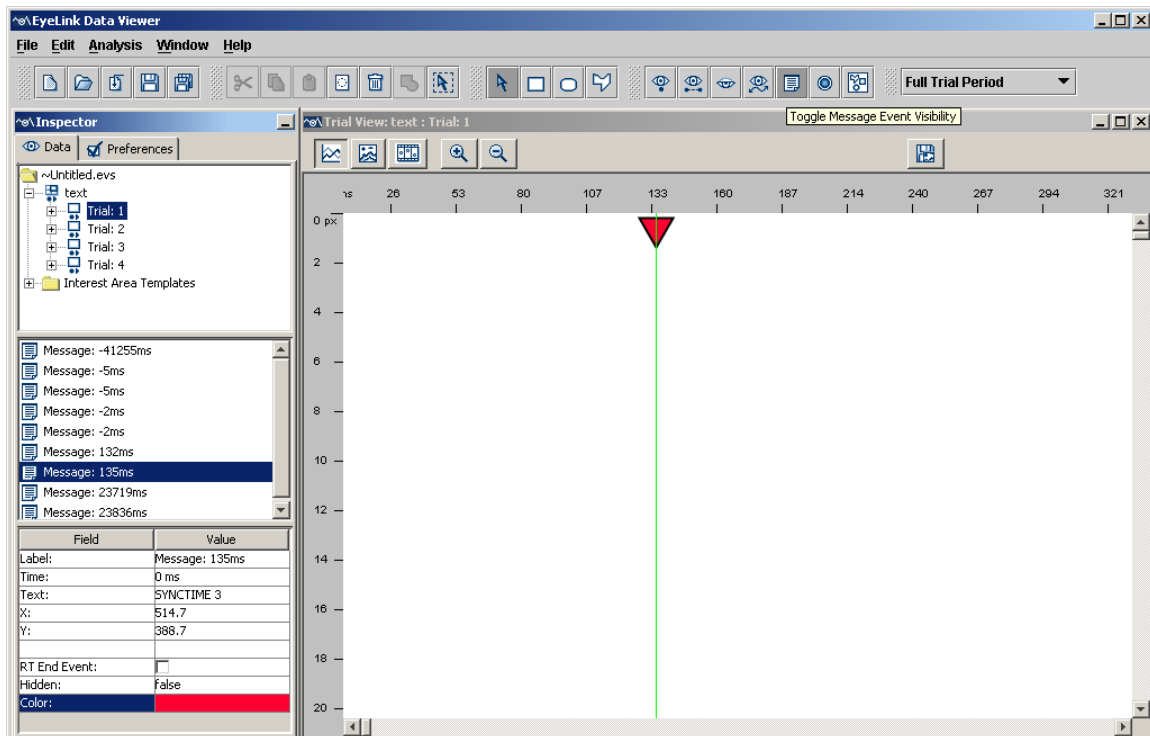


Figure 5-7. Sample Temporal Graph View of Message Events (Note that the default size of the message symbol in the temporal-graph preference settings has been changed).

5.5.1 Import Message List

Experiments should place certain messages into the EDF file to enable Data Viewer to process the files more efficiently. Examples of these commands include defining the image to overlay, specifying trial variables, defining a start time of RT definition, etc. These messages will be timestamped with an accuracy of 1 millisecond from the time sent, and can be used to mark important events such as display changes. If the user forgot to write these messages when the experiment was created, these messages can still be imported from a text file.

The loaded message list file must be a plain tab-delimited text file. Each line must be a valid EyeLink MSG line in the format of:

MSG EDF_time [Offset] Message_text

where:

- EDF_time must be an integer. If this field is positive, it is assumed to be the absolute EDF timestamp to be used for the message (e.g., a value of 62798072 for EDF_time field will place the message in a trial which started after 62798072 and ended before 62798072). If EDF_time is negative, the -EDF_time = trial index of the message in the EDF file. A value of -1 would mean the first trial in the file and -4 would mean the 4th trial in the file. In this case, the message EDF time is set to the start of the trial time in this case.


- Offset is optional, and must be an integer if specified. Offset is subtracted from EDF_time to generate the real message time. For example, a message line of “MSG 2435806 -14 !V APLAYSTART 0 1 waves/1.wav” suggests that the event the message was referring to (APLAYSTART) actual happened at time 2435820.
- Message_text can be any string up to 100 characters in size that starts with a non-numeric character.
- Lines that do not follow these rules are ignored.

e.g.,

```
MSG 2435806 -14 !V APLAYSTART 0 1 waves/1.wav
MSG 2436376 -14 !V APLAYSTOP 12578 1 waves/1.wav
MSG 62798072 0 POSN 174 384
MSG 62798072 0 POSN 185 384
```


To load the message list file:

- 1) Select the EDF file label in the treeview panel of the Inspector Window.
- 2) Click on the right mouse button and select “Import Message List”.
- 3) In the following “Open” dialog box, go to the directory where the message list file is stored and choose the file.

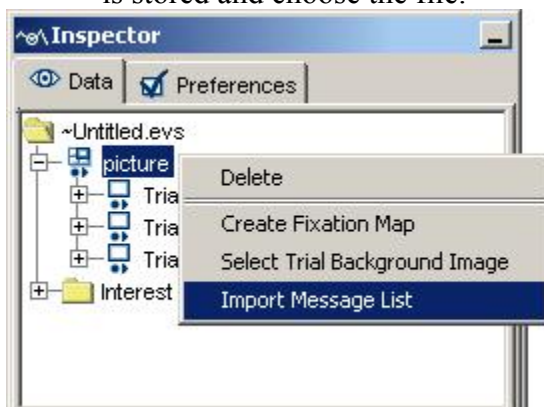

After the message list file is imported into the session, a dialog box will be presented to show the number of valid messages imported.

5.6 Button Events

Button events can be toggled on/off by clicking the Toggle Button Event Visibility button . A button event is represented as a colored triangle in the temporal graph view and a colored square in the spatial overlay view, centering on the current gaze position. The following table lists the properties of a button event. Those fields that are editable in the viewer are marked with an asterisk.

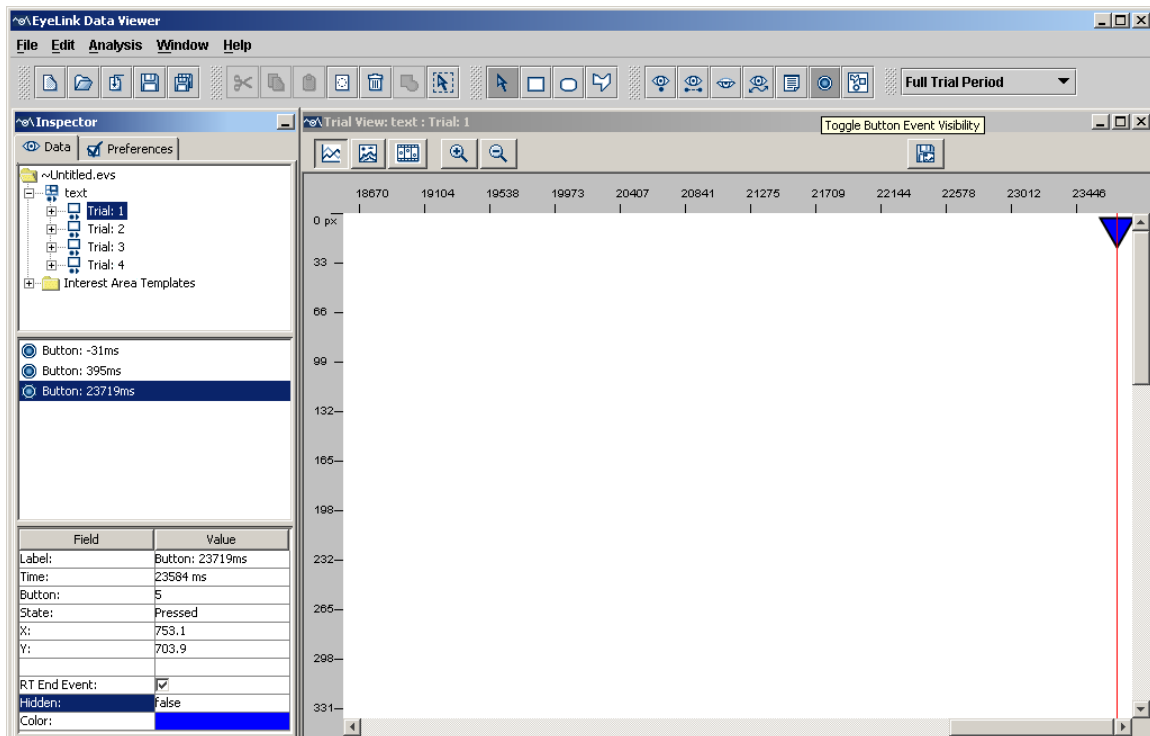



Figure 5-8. Sample Temporal Graph View Showing the Button Events (Note that the default size of the button symbol in the temporal-graph preference settings has been changed).

Field	Contents
Label *	Label of the button event
Time	Trial time when the button was pressed/released. Note that if the button event occurs before the trial starts, a negative time value will be displayed
Button	Button ID
State	Button pressed or released
X, Y	Concurrent gaze coordinates when the button is pressed/released (-100 if no gaze information was available for the time of the button event).
RT End Event *	Whether the current button event is the reaction time end event (see reaction time definitions).
Hidden	Whether the current button event is hidden from viewing and output report. This feature was disabled since version 1.6.
Color *	To achieve better visibility, the color of the currently selected button event can be changed by clicking on the color toolbox.

5.7 Samples

Samples can be toggled on/off by clicking the Toggle Eye Sample Visibility button  in the desktop toolbar. In the spatial overlay view (Figure 5-9, left panel), the samples are overlaid on the trial bitmap, forming the sample path of the viewing process. In the temporal graph view (Figure 5-9, right panel), the *xy* coordinates of the samples as well as the concurrent pupil size are plotted as a function of trial time.

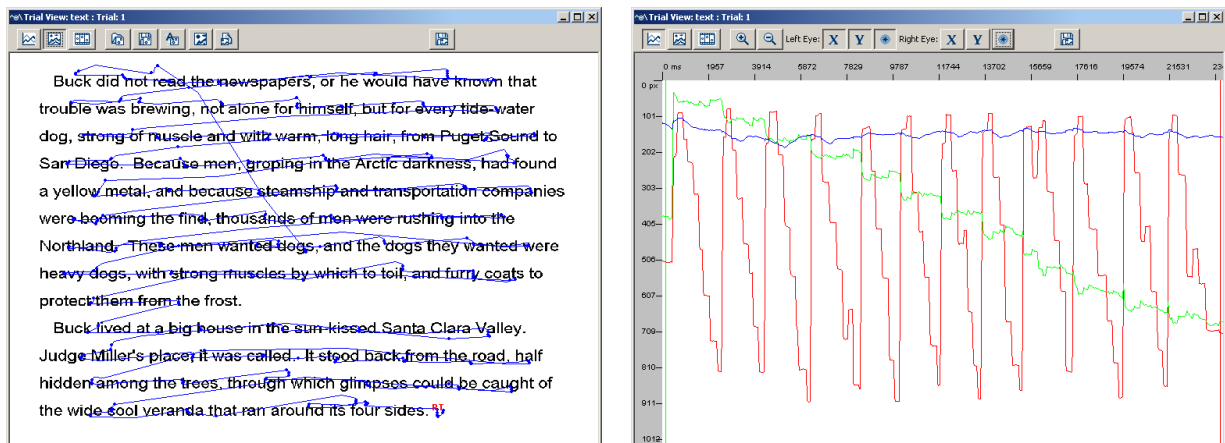



Figure 5-9. Sample Spatial Overlay View (left) and Temporal Graph View (right) of Raw Sample Data

Note: By default, samples are not loaded into the viewer when a data file is imported. To enable loading samples into the data viewer, change the Load Samples property in the Data Loading preference settings. If you want this change to be persistent over future viewing sessions, you may consider saving this as default setting. This can be done by selecting the topmost preference tree node, clicking the right mouse buttons and selecting "Save Properties as Defaults" (see Preference Settings).

Tip: In the Spatial Overlay trial view, if samples are visible at the same time as events, when an event is selected, all samples used to generate that event are also selected.

5.8 Interest Areas

To work on interest areas in a viewing session, make sure the **main** "Toggle Interest Area Visibility" button () in the application toolbar is turned on. To access interest area data for individual trials, first expand the trial by clicking on the + next to it in the tree view Panel of the Inspector window, then click on the Toggle interest area visibility button for that trial. The node will be labeled "Custom Interest Area Set" if custom interest areas are loaded or created or "Empty Interest Area Set" if no interest areas are defined. The middle panel of the inspector window lists all of the available interest areas

for the currently selected trial. The lower panel shows the properties of the selected interest area.

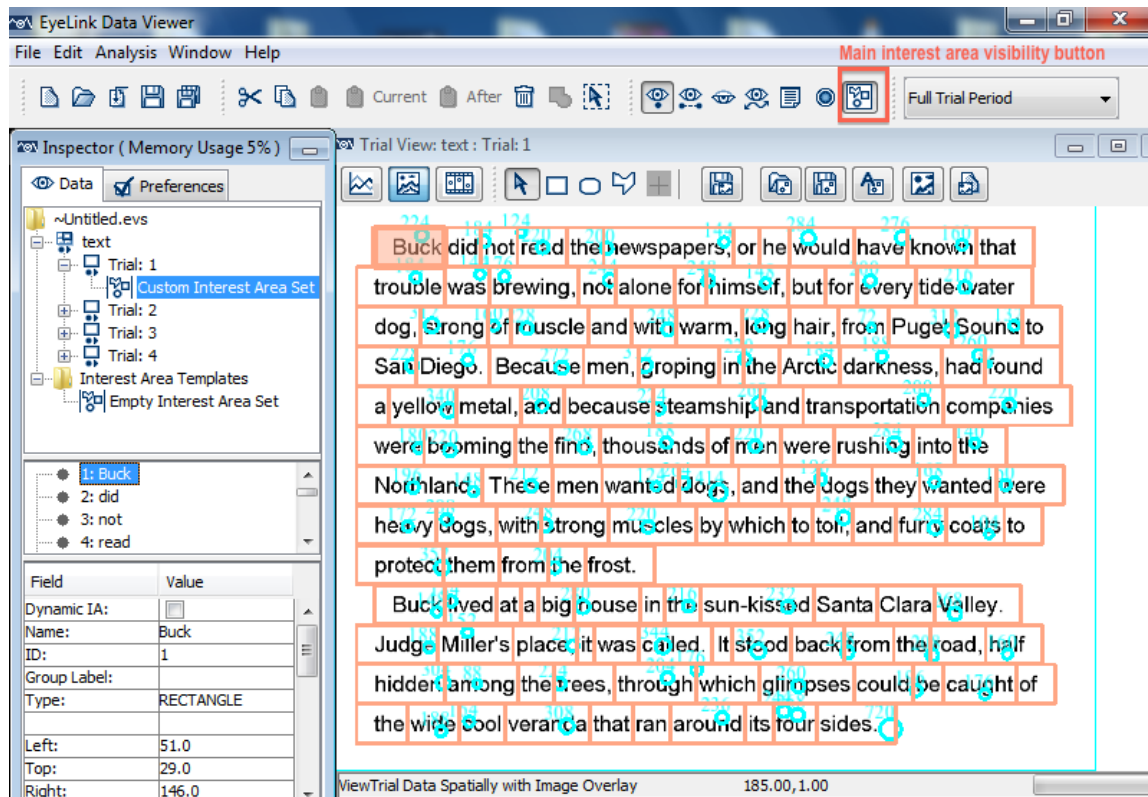


Figure 5-10. Sample Text Display Overlaid with Interest Areas

5.8.1 Format of static and dynamic interest areas

Version 2.0 of Data Viewer allows both static and dynamic interest areas. Static interest areas either have no time stamp (in which case they are present for the entire duration of the trial or selected Interest Period) or a single time stamp which indicates the time of their onset relative to the beginning of the trial. Dynamic interest areas have two time stamps - the first indicating the time of their onset, and the second indicating the time of their offset.

Thus interest areas can be encoded in one of the following three formats:

Type 1:

RECTANGLE <id> <left> <top> <right> <bottom> [label]

ELLIPSE <id> <left> <top> <right> <bottom> [label]

FREEHAND <id> <x1, y1> <x2, y2> ... <xn, yn> [label]

e.g., RECTANGLE 1 0 0 512 384 images\hearts.jpg

Type 2:

```

start_offset RECTANGLE <id> <left> <top> <right> <bottom> [label]
start_offset ELLIPSE <id> <left> <top> <right> <bottom> [label]
start_offset FREEHAND <id> <x1, y1> <x2, y2> ... <xn, yn> [label]
e.g., -500 RECTANGLE 1 0 0 512 384 images\hearts.jpg

```

Type 3:

```

start_offset end_offset RECTANGLE <id> <left> <top> <right> <bottom> [label]
start_offset end_offset ELLIPSE <id> <left> <top> <right> <bottom> [label]
start_offset end_offset FREEHAND <id> <x1, y1> <x2, y2> ... <xn, yn> [label]
e.g., -500 -1500 RECTANGLE 1 0 0 512 384 images\hearts.jpg

```

The first two examples are static interest areas whereas the third example is a dynamic interest area. The "start_offset" and "end_offset" values are usually negative, indicating a future time relative to the start of the currently selected interest period (the default Full Trial Period will be used if no interest periods have been created for the viewing session). For example, if the current interest period starts at 800 ms following the trial start, the particular interest area of the above Type 3 example would start at 1300 ms and end at 2300 ms following the trial start. If the interest areas are read from a file that a message in the EDF file specifies (e.g., the "!V IAREA FILE" message in the EDF file), the "start_offset" and "end_offset" are relative to the time stamp of that message in the EDF file.

As with other events in a viewing session (such as fixations and saccades), interest areas are displayed in the event list. Static interest areas are represented by a dot and dynamic ones are represented by a folder; there will be a separate folder for each of the dynamic interest areas in a trial. Each dynamic interest area may contain several instances to describe the position of the interest area across different time segments (a dynamic interest area object must contain at least one instance). For example, a fixation screen (with a "FIXATION_SCREEN" message written to the EDF file) is shown for 1000 ms before a target screen (with a "SYNCTIME" message) is presented with a red object appearing at (674, 546, 774, 646). The object moves to (462, 634, 562, 734) at time 800 ms after its onset, to (249, 546, 349, 646) at 1600 ms, and to (162, 334, 262, 434) at time 2400 ms and stays there for 800 ms before a blank screen is shown (with a "BLANK_SCREEN" message). A dynamic interest area of four instances can be coded (shown below) to describe the behavior of the red object, assuming that an interest period is created starting from the SYNCTIME message and this interest period is selected.

0	-799	RECTANGLE2	674	546	774	646	red
-800	-1599	RECTANGLE2	462	634	562	734	red
-1600	-2399	RECTANGLE2	249	546	349	646	red
-2400	-3199	RECTANGLE2	162	334	262	434	red

You will see the following four entries in the "2:red" folder of the event list, representing four instances of the dynamic interest area (2:red). See section 24 for a summary of the various methods available for creating dynamic interest areas in a viewing session.

2:red (0:799)

2:red (800:1599)
 2:red (1600:2399)
 2:red (2400:3199)

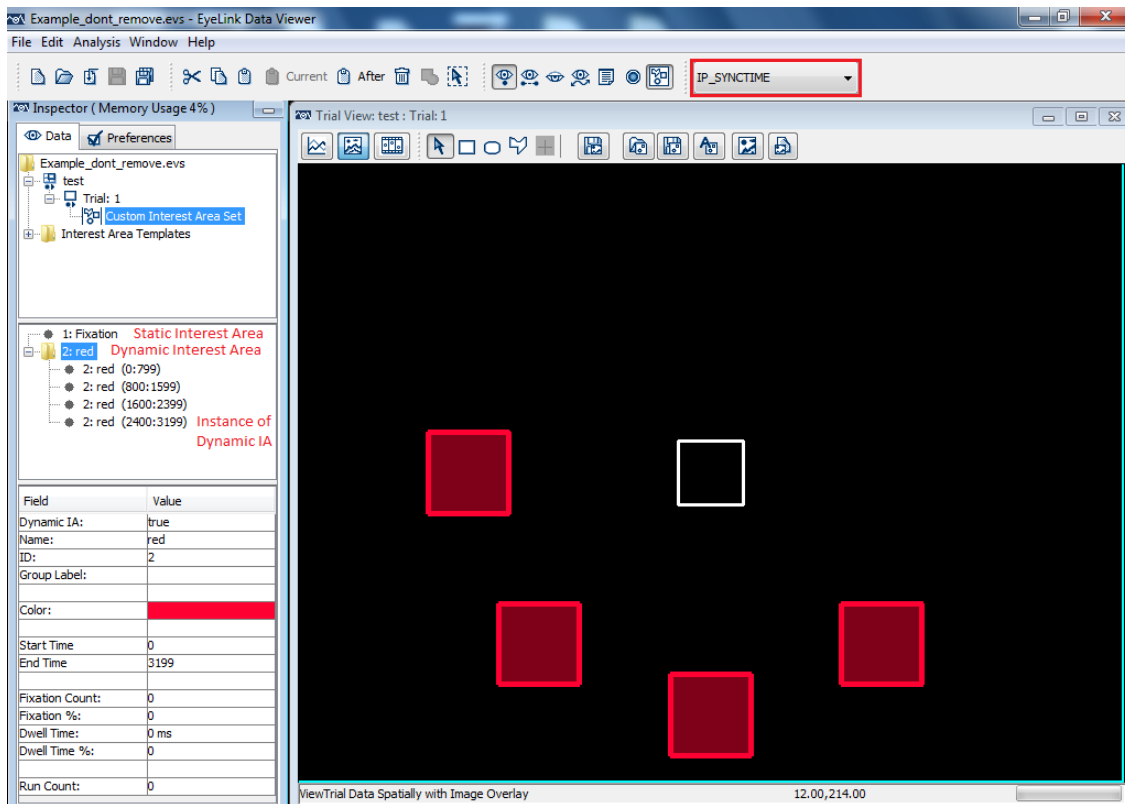


Figure 5-11. Sample Interest Area Instances in One Interest Period

The start time and end time of the instances will be different if you then change the interest period. For example, the start time and end times will look like the following if you create and select a new interest period starting from the “FIXATION_SCREEN” message time.

-1000	-1799	RECTANGLE2	674	546	774	646	red
-1800	-2599	RECTANGLE2	462	634	562	734	red
-2600	-3399	RECTANGLE2	249	546	349	646	red
-3400	-4199	RECTANGLE2	162	334	262	434	red

You will see the following four entries in the “2:red” folder of the event list, representing four instances of the dynamic interest area (2:red) under the new interest period.

2:red (1000:1799)
 2:red (1800:2599)
 2:red (2600:3399)
 2:red (3400:4199)

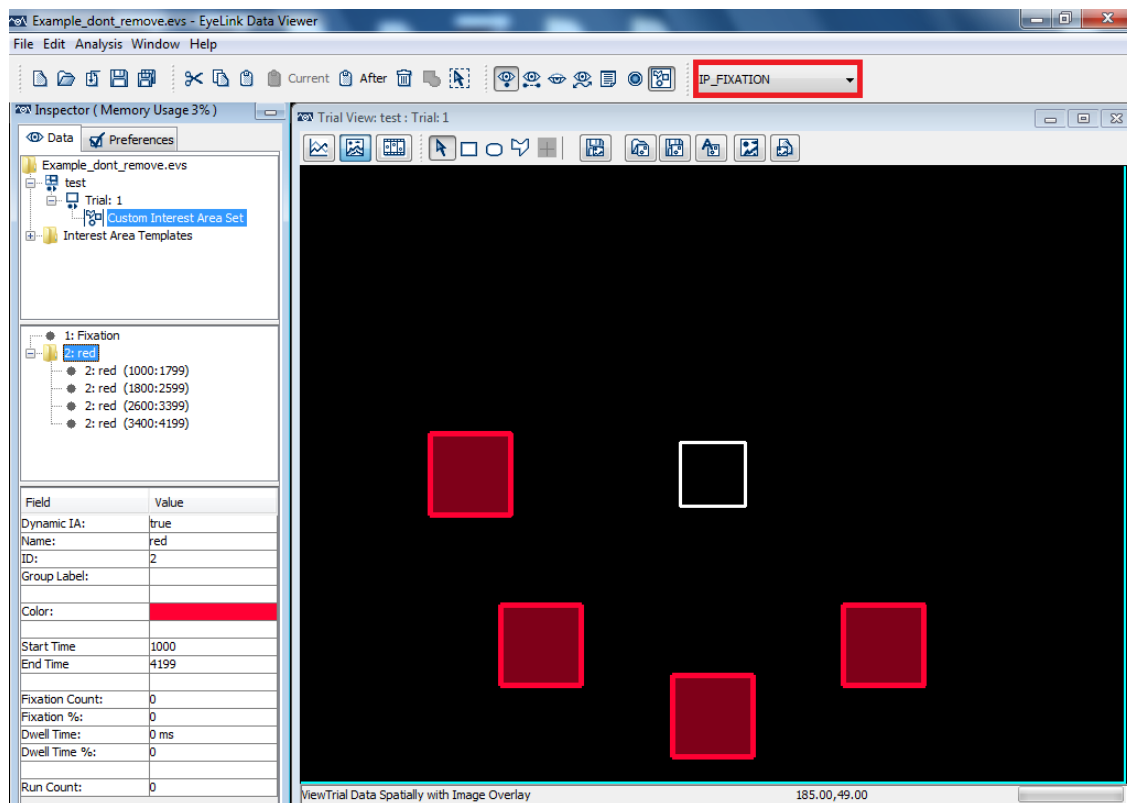


Figure 5-12. Sample Interest Area Instances in Another Interest Period

Clicking on the instance in the event list will highlight the interest area in the spatial overlay view and vice versa. If the trial view window is set to the animation view, clicking on the instance will set the playback time of the animation view to the start time of the instance if the target interest area is not currently displayed.

5.8.2 Interest Area Properties

The following table lists the properties of an interest area. Those fields that are editable in the viewer are marked with an asterisk (*).

Property	Property Description
Dynamic IA	Whether the current interest area is a dynamic one (checked) or a static one (unchecked). This field is not directly editable. You can convert a static interest area into a dynamic one or vice versa through the “Convert to DynIA/StaticIA” option in the popup menu when you right click on the interest area.
Name *	Text of the interest area.
ID	Ordinal ID of the interest area.
Group	Label of the group to which the current interest area belongs.
Type	Type of interest area (rectangular, elliptical, or freehand).




Left *	X-coordinate of the left-most edge of a rectangular or elliptical interest area.
Top *	Y-coordinate of the top-most edge of a rectangular or elliptical interest area.
Right *	X-coordinate of the right-most edge of a rectangular or elliptical interest area.
Bottom *	Y-coordinate of the bottom-most edge of a rectangular or elliptical interest area.
Center	Pixel coordinate (x, y) for the center of the interest area.
Pixel Area	Area of the interest area in pixels.
Color *	To achieve better visibility, the color of the interest area or its instances can be changed through a color selector.
Time	EDF file time of the message specifying the interest area if it was created using the '!V IAREA' command during recording. If the interest area was created manually or from an interest area template file during the viewing session, the time value will be -1 and this attribute will not be visible from the property table. This time stamp is useful in case you need to filter interest areas based on the currently-selected interest period (see “Show Interest Areas Pre Interest Period” and “Show Interest Areas Post Interest Period” preferences in “Preferences -> Data Filters”). This field is for static interest areas only.
Start Time relative to IP / End Time relative to IP	The start time / end time (in milliseconds) of the dynamic interest area or its instance relative to the start of the currently selected interest period. This field is for dynamic interest areas only. After instances of dynamic interest areas are created in the viewing session or imported from an interest area file, the start time and end time will be adjusted if a new interest period is chosen.
Start Time relative to Trial / End Time relative to Trial	The start time / end time (in milliseconds) of the dynamic interest area or its instance relative to the start of the trial. This field is for dynamic interest areas only and its value remains fixed regardless of the interest period that is currently selected.
Fixation Count	Total number of fixations within the interest area in the current interest period.
Fixation %	Percentage of total fixations in a trial falling within the current interest area in the current interest period.
Dwell Time	Total time (in milliseconds) spent on the current interest area in the current interest period.
Dwell Time %	Percentage of trial dwell time spent on the current interest area in the current interest period.
Run Count	Number of times the interest area was entered and left in the current interest period.

Note: The coordinate information (Left, Top, Right, Bottom, and Center) is applicable to rectangular and elliptical interest areas only.


5.8.3 Manually Creating Interest Areas

To create an interest area:


- 1) Select the trial view window by clicking on or in it, so that the window is active.
- 2) Select the Interest Area shape that you want to create from the toolbar of the trial view window:

	Rectangular Interest Area Shape is active
	Ellipse Interest Area Shape is active
	FreeHand Interest Area Shape is active

For rectangle and ellipse interest areas:

- 3) Hold down the left mouse button in the top-left corner of where you want the interest area to be created.
- 4) Drag the mouse down to the right until the selection region covers the area that you want to select.
- 5) Release the mouse button.
- 6) In the “New Area of Interest Data” dialog box, fill out the “label” field. If you are creating a static interest area, leave the “Dynamic IA” box unchecked and press Enter. For a dynamic interest area, you will need to further configure its properties - please see the subsection on configuring dynamic interest areas for details.
- 7) Repeat steps 3-6 to create more interest areas of the same type. Press the Selection mode icon () from the toolbar of the trial view window to stop creating interest areas.

For FreeHand interest area shapes:

- 3) Click the left mouse button on the first point of the shape you want to create.
- 4) Move the mouse to the second point and click the left mouse button again. A line joining the points will be displayed.
- 5) Repeat step 4 for each point of the freehand shape.
- 6) Press the ENTER key when you have specified all points you wish to create. Data Viewer will automatically join the first and last points of the interest area. Note that you can cancel the creation of a freehand interest area by pressing the ESC key.
- 7) In the “New Area of Interest Data” dialog box, fill out the “label” field. If you are creating a static interest area, leave the “Dynamic IA” box unchecked and press Enter. For a dynamic interest area, you will need to further configure its properties - please see the following section for details.
- 8) Repeat steps 3-7 to create more interest areas of the same type. Press the Selection mode icon () from the toolbar of the trial view window to stop the creation process.

Note: Please avoid creating self-intersecting freehand interest areas, as this will cause problems for the proper calculation of pixel area.

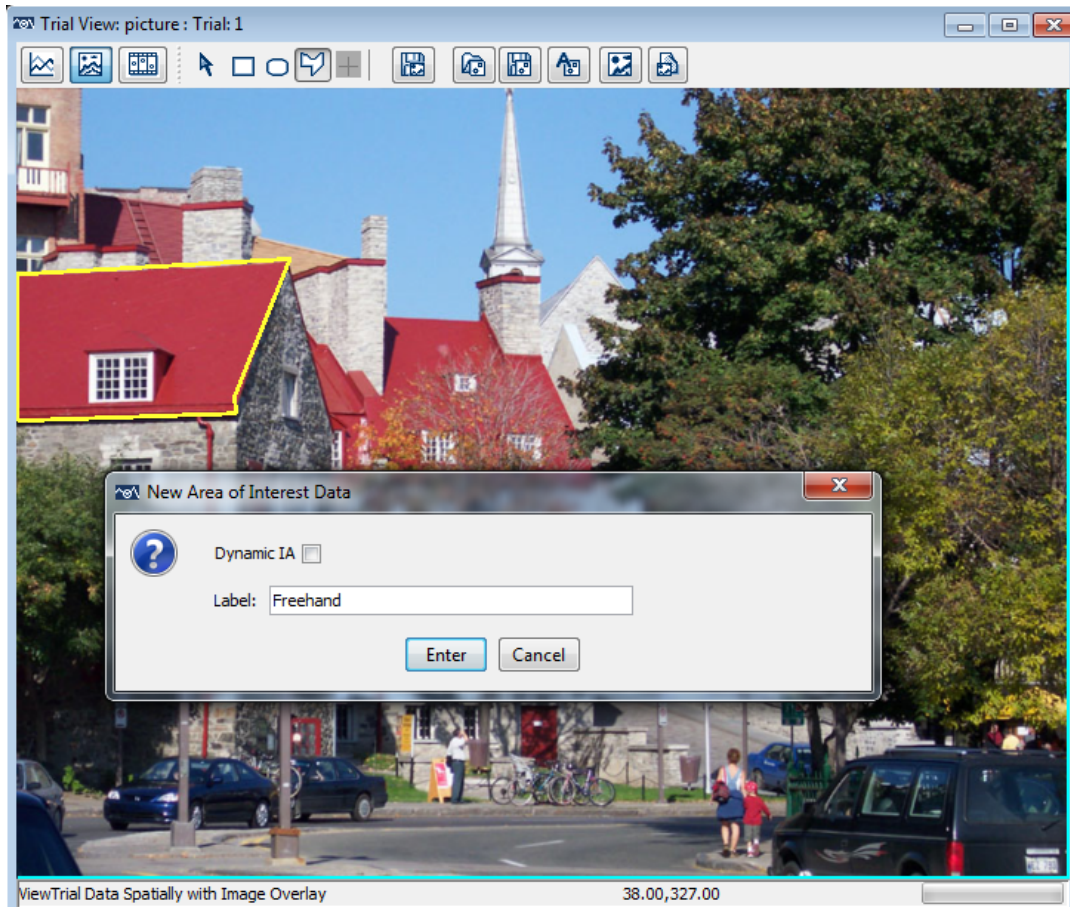


Figure 5-13. Creating a Freehand Interest Area

Configuring Dynamic Interest Areas:

If the “Dynamic IA” checkbox is ticked additional properties will be shown in the New Area of Interest Data dialog:

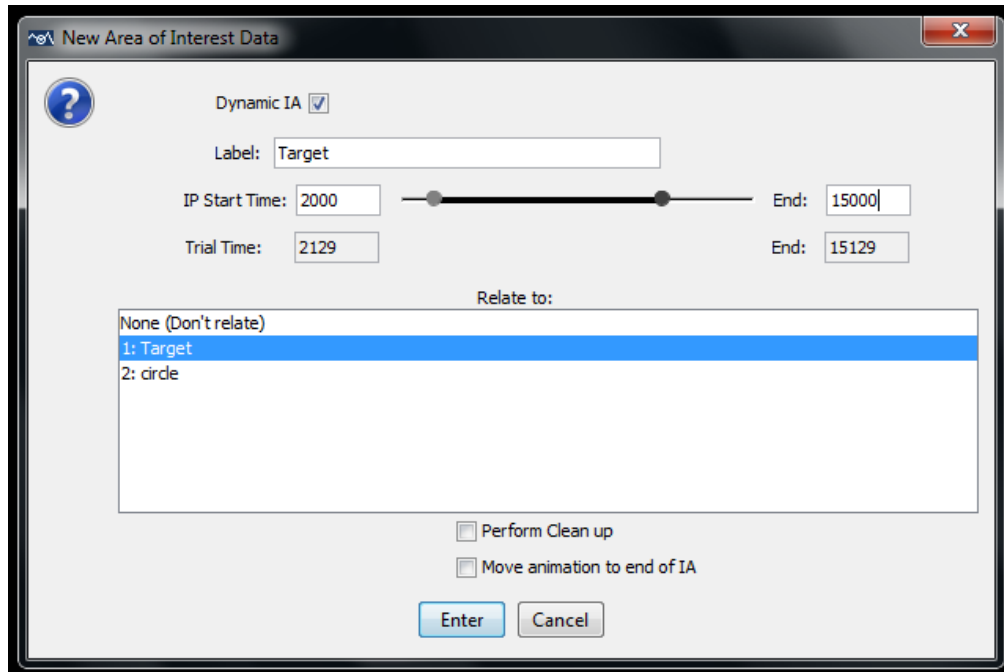



Figure 5-14. Sample Interest Area Instances in Another Interest Period

- 1) IP Start Time: The start time (in milliseconds) of the current instance of the interest area relative to the start of the currently selected interest period. **It is critical that an interest period has been created and selected before creating a dynamic interest area.** You can either type in the value in the edit box or click on the left indicator on the sliding bar and drag it to the intended position.
- 2) IP End Time: The end time (in milliseconds) of the current instance of the interest area relative to the start of the interest period. You can either type in a value in the edit box or click on the right indicator on the sliding bar and drag it to the intended position.
- 3) Trial Time: The start time (in milliseconds) of the current instance of the dynamic interest area relative to the start of the trial. This field is read-only and changes along with the “IP start time”.
- 4) Trial End Time: The end time (in milliseconds) of the current instance of the dynamic interest area relative to the start of the trial. This field is read-only and changes along with the “IP end time”.
- 5) “Relate to”: this specifies whether the current instance is a member of a new interest area (“None (Don't Relate)”) or a new instance of an existing dynamic interest area. If the latter is the case, select the existing interest area from the list below. The newly-created instance will be created in the folder of the selected interest area in the event list.
- 6) Perform Clean up: If this option is checked, Data Viewer will perform a Clean up operation on the instances of the interest area (e.g., to trim instances with overlapping times - see section “13 Cleaning Up Dynamic Interest areas” below).
- 7) Move animation to end of IA: If this option is checked, the playback time of the

animation view will be set to 1 millisecond after the end point of the newly created interest area instance.


5.8.4 Creating Interest Area: Image Segmentation

If a trial bitmap has already been loaded, the Data Viewer can create interest areas for the trial by auto-segmentation. This is done by clicking the “Auto Segment Image to Interest Area” button () in the toolbar of the spatial overlay view window.

Note: Using the auto-segmentation feature will remove all of the existing interest areas for the trial. Auto segmentation is more feasible for those displays in which the individual items can be easily isolated, such as in reading and visual search studies where a background of a single color is present, but not good for some visual scenes in which the boundaries between individual elements are not clearly defined (see Section 8.7 on preference settings for image segmentation).

5.8.5 Saving Interest Areas


Interest areas that have been created for any one trial can be saved into a template file and re-used on other trials. To save interest areas:

- 1) Click the **Save the Interest Area Set to Disk** button () on the toolbar of the trial view window.
- 2) In the Save dialog box, go to the directory where the interest area set should be saved.
- 3) In the File name text box, type in the name of the file. Data Viewer will automatically fill in the file extension (.ias).
- 4) Press the Save button.

In the saved interest area file, static interest areas will not have a time stamp. For dynamic interest areas, the time stamps associated with each instance of the interest area will be based on the currently selected interest period. So it is critical that the interest period is set correctly before saving the template. Please note that the time stamps will change if an incorrect interest period is selected when saving – this will mean that when the template is applied to other trials (when the interest period is set properly) the instances will not have the correct time stamps.

5.8.6 Importing Interest Area File (for a specific trial)

Interest areas from a file can be imported and applied to a specific trial.

- 1) Click the **Import Interest Area File** button () on the toolbar of the trial view window
- 2) In the Load dialog box, browse to the location of the interest area file that you want to import and select it.
- 3) Click Load.

The interest areas imported from the file will be added on top of the existing interest areas in the trial. For dynamic interest areas, the start and end time will be mapped relative to the currently selected interest period. If no interest period is created for the session, the default “Full Trial Period” will be used.

5.8.7 Emptying an Interest Area Set

To empty the current interest area set for a trial:

- 1) Click on the target trial in the tree view whose interest areas are to be emptied
- 2) In the property panel of the trial, find the entry “Interest Area Set” and choose “Empty Interest Area”.

Existing interest areas in a recording session file or a group of trials can be similarly cleared by setting the “Default Interest Area Set” entry of the property table to “Empty Interest Area Set”.

5.8.8 Loading Interest Area Templates

If the same interest areas are used across a group of trials, instead of repeating the procedure of creating interest areas for each individual trial, an interest area template can be used to supply the interest area information to a group of trials. To import an interest area template file into a viewing session:

- 1) From the application menu, choose:
 File -> Import Data -> Interest Area Template
- 2) In the Load dialog box, browse to the location of the interest area (.ias) file(s) that you want to import and select it/them.
- 3) Click Load.

The interest area template files imported will be listed in the “Interest Area Template” section, which is at the bottom of the tree view in the Inspector window.

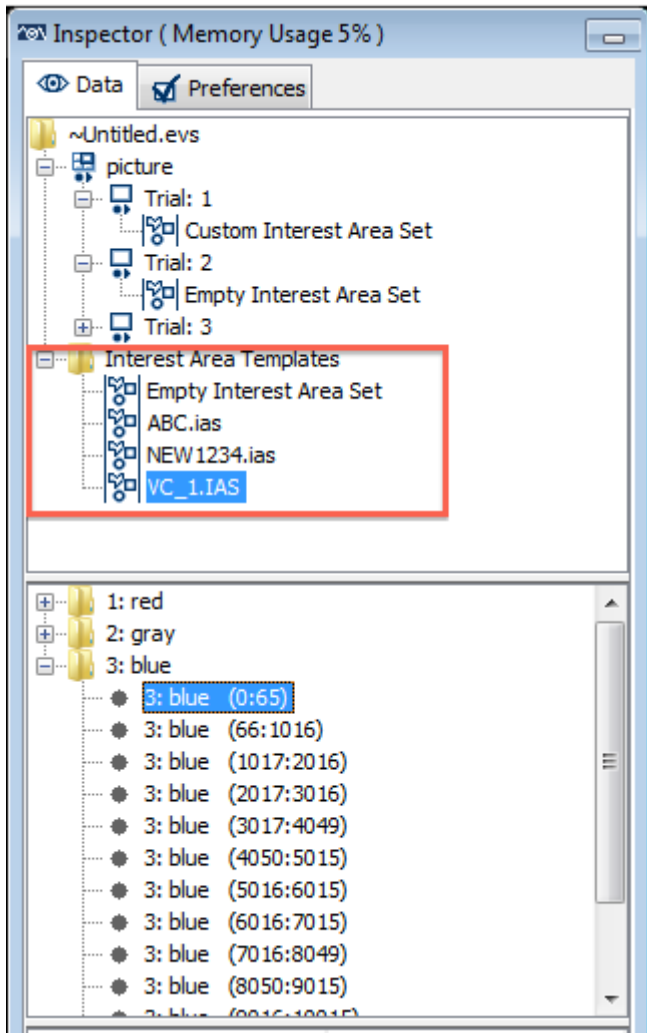


Figure 5-15. Interest Area Templates Listed in the Tree View

5.8.9 Using Interest Area Templates

An interest area template, if already loaded into the viewing session, can be used to apply the interest area data across a group of trials (see section 6.1.3 for information on the trial grouping feature of Data Viewer). To apply an interest area template to a trial or group of trials,

- 1) Click on the viewing session icon (📁) and perform trial grouping if necessary.
- 2) In the tree view panel of the data tab, select the trial/group onto which the interest area template should be applied.
- 3) In the property panel of the target trial or trial group, find the entry “Interest Area Set” and select the intended interest area template from the dropdown list.


Important: When applying an interest area template to a trial or a group of trials, the start time and end time of the dynamic interest areas that are applied to the trial/group of trials

will be set relative to the currently selected interest period. If no interest period is created for the viewing session, the default “Full Trial Period” will be used. So it is very important that a proper interest period be created and selected before applying interest area templates containing dynamic interest areas to a trial or group of trials!





5.8.10 Deleting Interest Area Templates

Interest area templates can be deleted from the template list. In the tree view panel of the Inspector window, find the "Interest Area Templates" folder where all of the interest area templates are located. Select the target interest area template from the list, click the right mouse button, and select "Delete". If the deleted interest area template is used by a trial, the "Interest Area Set" property of that trial will be set to "Empty Interest Area Set" and existing interest areas will be removed.

5.8.11 Selecting Interest Areas

To select an interest area or an instance of a dynamic interest area, place the mouse cursor over the interest area in the tree view window and click the left mouse button. The interest area is highlighted in both the trial view window and the event list. To select more than one interest area, hold down the CTRL key on Windows or the Command key ⌘ on Mac OS X and select the target area by left-clicking - repeat until all of the desired interest areas are selected. If all of the interest areas in a trial are to be selected, simply click the **Select All Visible Items** button () on the toolbar (or click the right mouse button and select “Select All”).

Selection can also be made from the event list in the inspector window by clicking on the “Custom Interest Area Set” node of the trial in the tree view and selecting the interest areas from the event list panel. The selected interest area will be highlighted in the trial view window. When an instance of a dynamic interest area is selected from the event list, the playtime of the animation view will be set to the start time of that instance.

Note: If you are not able to select any interest area or instance, please make sure none of the interest area shape buttons (, , or ) are active. You may press the Selection mode icon () from the toolbar of the trial view window to stop creating interest areas.

If there are eye events within the interest area to be selected, clicking on the interest area may end up selecting the eye events instead. To make a more precise interest area selection, you may toggle off the visibility of the eye events first, make the interest area selection, and then toggle the visibility of the eye events back on if desired.

5.8.12 Copying and Pasting Interest Areas

A new interest area can be created by copying and then pasting an existing one. Select the target interest area, click on the right mouse button, choose the “Copy” option from the

popup menu, and then choose the “Paste” option. This procedure can be used to create a copy of an existing static interest area, the entire dynamic interest area folder, or an instance of a dynamic interest area.

You will see additional “Paste Current” and “Paste After” option after copying an instance of a dynamic interest area. The “Paste Current” option, applicable to the animation view only, will create a new interest area instance with the start time being the current playtime of the animation view. The duration of the new instance is the same as the original instance. The “Paste After” option is available in both the animation view and the spatial overlay view. This option will create a new interest area instance of the same duration immediately following the end time of the original one.

5.8.13 Cleaning Up Dynamic Interest Areas


When creating dynamic interest areas, you may occasionally create instances of an interest area that overlap in time. The overlapping interest area instances can pose a problem for data analysis (e.g., when calculating dwell time on a dynamic interest area). Data Viewer provides a clean up option so that you can fix any overlapping timing issues for an interest area, a trial, a group of trials, or the entire viewing session. When cleaning up two interest area instances with overlapping times, if the two instances start at the same time, the more recently-created one will be preserved; otherwise, the overlapping portion of the instance that starts earlier in time will be removed. For example, instance A (start time a1 and end time a2) starts before instance B (start time b1 and end time b2). The cleanup process will preserve the original instance B while trimming the original instance A so that only the portion from a1 to b1-1 remains.



5.8.14 Mouse Record Mode

Mouse Record mode provides an easy interface to create dynamic interest areas in the animation view. This mode allows you to first make a copy of an existing interest area instance, and then insert subsequent new instances with a mouse click at the intended playback time.

- 1) Select an instance of a dynamic interest area either in the tree view or by clicking on it in the trial view window.
- 2) Click the right mouse button, and choose the “Mouse Record” mode option from the popup menu (see the screen capture below).
- 3) Now set the trial view window to the animation view
- 4) Use the slider bar at the top of the trial view window to set the animation view to the intended playback time (you may use the left and right cursor key to step backward and forward in playback time).

- 5) Now move the mouse within the trial view window (note that the mouse cursor changes into a +). Click the left mouse button on top of the intended location – a new instance will be created following the mouse click.
- 6) Repeat this step until all of the instances are created.
- 7) End the Mouse record mode by clicking on the Selection mode icon () from the toolbar of the trial view window, or by pressing the ESC key.

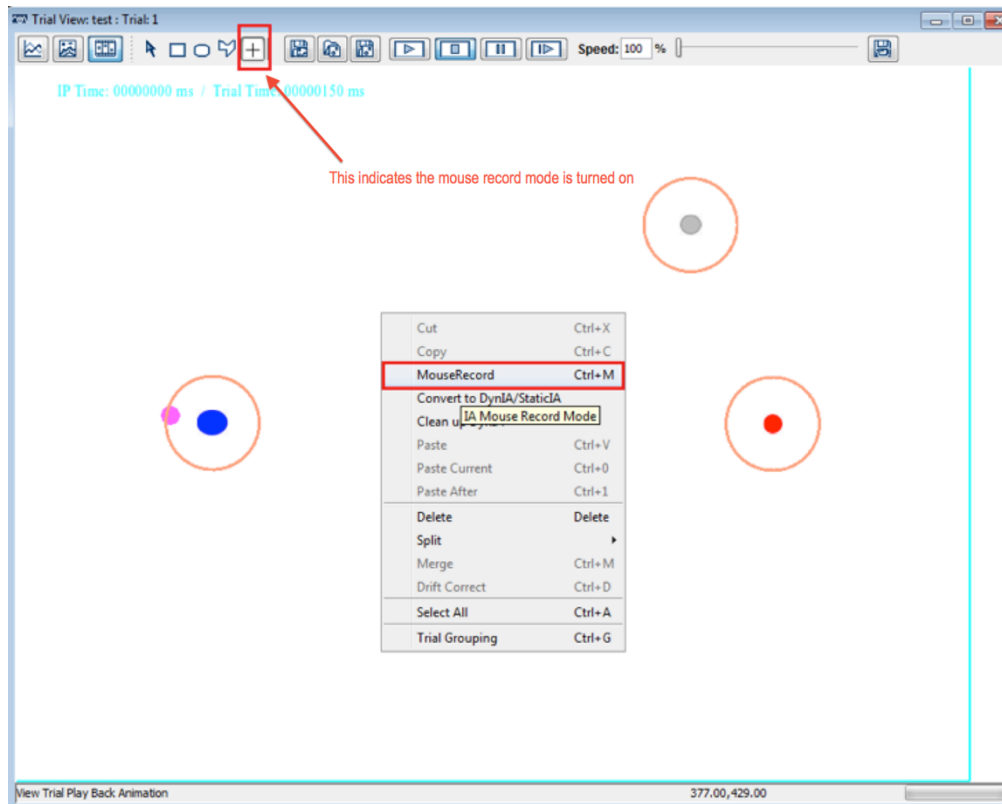


Figure 5-16. Mouse Record Mode

Note: although you can create interest areas while the animation is playing back in the animation view, it is probably a better practice to set the exact time position and create the interest areas at that specific time point. Creating the interest areas while the playback is in progress can result in lags in the start time of the interest areas.

5.8.15 Moving an Interest Area

Selected interest area(s) can be moved by following the steps below:


- 1) Select the interest area(s) (make sure the selection is highlighted in the trial view window).
- 2) Hold down the ALT key.
- 3) Place the mouse cursor on top of the interest area(s) and press and hold the left mouse button.
- 4) Drag the interest area(s) to the intended location.

- 5) Release the mouse button.

The position of the selected interest area(s) can also be adjusted by pressing the left, right, up, and down keys while holding down the ALT key.

5.8.16 Resizing Rectangular and Elliptical Interest Areas

A rectangular or elliptical interest area can be resized by following these steps:

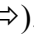- 1) Select the interest area (make sure the selection is highlighted in the trial view window).
- 2) Hold down the ALT key.
- 3) Place the mouse over the edge to be dragged (the arrow cursor changes into a sizing cursor ).
- 4) Press the left mouse button and drag the edge to the desired location.
- 5) Release the mouse button.


To carry out a more precise adjustment for a rectangular or elliptical interest area:

- 1) Select the interest area to move or adjust.
- 2) In the property panel of the interest area, select the entry for “Left” and type in the desired x-coordinate of left-most edge.
- 3) Select the entry for “Top” and type in the desired y-coordinate of top-most edge.
- 4) Select the entry for “Right” and type in the desired x-coordinate of right-most edge.
- 5) Select the entry for “Bottom” and type in the desired y-coordinate of bottom-most edge.

Please note that the coordinates must be made such that the “Left” value will be smaller than the “Right” value and the “Top” value will be smaller than the “Bottom” value.

5.8.17 Resizing Freehand Interest Areas


A freehand interest area can be resized by following these steps:

- 1) Select the interest area (make sure the selection is highlighted in the trial view window).
- 2) Hold the ALT key down.
- 3) Place the mouse over the vertex to be moved (the move cursor changes into a hand cursor  and a red circle is drawn on top of the vertex).
- 4) Press the left mouse button and drag the vertex to the desired location.
- 1) Release the mouse button.

5.8.18 Adding a Point to a Freehand Interest Area

To add a new point to a freehand interest area:

- 1) Select the interest area (make sure the selection is highlighted in the trial view window)

- 2) Hold the ALT key down.
- 3) Place the mouse over the existing edge where a point will be added (the arrow cursor changes into a hand cursor )
- 4) Press the left mouse button. A new vertex will be created.
- 5) Now move the newly added vertex to the intended location.

5.8.19 Deleting an Existing Point from a Freehand Interest Area



To remove an existing point from a freehand interest area:

- 1) Select the interest area (make sure the selection is highlighted in the trial view window)
- 2) Hold the ALT key down.
- 3) Place the mouse over the vertex to be removed (the arrow cursor changes into a hand cursor and a red circle is drawn on the vertex).
- 4) Click on the right mouse button, and choose the “Delete” option from the popup menu.
- 5) When the delete confirmation dialog box is displayed (“Do you want to delete the selected point from the Freehand Interest Area”), choose “Yes”. Data Viewer will automatically join the two points around the deleted one.

5.8.20 Deleting Interest Areas

To delete interest areas:

- 1) First toggle off all of the eye event visibility buttons in the application toolbar so that interest areas can be deleted without removing any eye events within them.
- 2) Select the interest areas to be deleted.
- 3) Press the right mouse button.
- 4) Select the Delete option from the popup menu.
- 5) Confirm that the IAs should be deleted.

Interest areas can also be deleted by clicking the Delete Selected Items button () or the Cut Selected Item button () on the application toolbar.

5.8.21 Splitting Interest Areas

Rectangular and elliptical interest areas can be split either horizontally or vertically. To split an interest area:

- 1) Select the interest area to split.
- 2) Press the right mouse button.
- 3) Select “Split” and then “Horizontally” or “Vertically” from the popup menu.

When splitting an instance of a dynamic interest area, the duration of the instance is split as well - one of the newly created instances takes the first half of the original duration and the other takes the second half of the duration.

5.8.22 Merging Interest Areas / Instances of a Dynamic Interest Area

Two selected rectangular or elliptical interest areas can be merged in the following way:

- 1) Click "Toggle Fixation Event Visibility" button (👁️) on the application toolbar to turn off the fixation visibility so that fixation events will not be merged.
- 2) Press the CTRL key on Windows or the Command key ⌘ on Mac OS X to select the interest areas to be merged.
- 3) Click the right mouse button to bring up a popup menu, select "Merge" from the list.

Multiple instances of a dynamic interest area can also be merged. The start time of the newly created instance will be the start of the earliest instance and the end time will be end of the latest instance

5.8.23 Conversion between a Static Interest Area and a Dynamic Interest Area


Unlike a static interest area, a dynamic interest area consists of one or multiple instances, each of which has a start time and end time. Data Viewer supports the conversion between these two types of interest areas: instances of a dynamic interest area can be converted into static interest areas by removing the start time and end time whereas a static interest area can be converted into a dynamic one, with the start time being 0 and end time being the duration of the trial. The conversion process preserves the spatial coordinate of the original interest area.

5.8.24 Examples of Creating Dynamic Interest Areas in a Viewing Session

This section briefly summarizes various methods of creating dynamic interest areas using the example given at the beginning of the Chapter. A fixation screen (with a "FIXATION_SCREEN" message written to the EDF file) is shown for 1000 ms before a display screen (with a "SYNCTIME" message) is presented with a red object appearing at (674, 546, 774, 646). The object moves to (462, 634, 562, 734) at time 800 ms after its onset, to (249, 546, 349, 646) at 1600 ms, and to (162, 334, 262, 434) at time 2400 ms and stays there for 800 ms before a blank screen is shown (with a "BLANK_SCREEN" message). A dynamic interest area of four instances can be coded (shown below) to describe the behavior of the red object, assuming that we create an interest period starting from the onset of the SYNCTIME screen.

0	-799	RECTANGLE2	674	546	774	646	red
-800	-1599	RECTANGLE2	462	634	562	734	red
-1600	-2399	RECTANGLE2	249	546	349	646	red
-2400	-3199	RECTANGLE2	162	334	262	434	red

There are several possible ways to create the instances of this dynamic interest area. Regardless of the method chosen, the first step would be creating an interest period and selecting it as the active interest period.

- 1) Click the "Interest Period List" combo box at the far-right hand of the application menu bar and select "Edit ...".
- 2) Click the *New* button () on the interest period manager dialog box.
- 3) In the general tab, enter a label for the current interest period in the label edit box and choose "Message Event" in the "Type" drop-down list.
- 4) Now click on the "Details" tab. Set the "Start Time Message Text" as "SYNCTIME" and "End Time Message Text" as "BLANK_SCREEN" (both message texts are case-sensitive without quotation marks).
- 5) Close the Interest Period Editor dialog. The "Interest Period List" combo box should now display the label of the newly created interest period.

Next, create the first instance of the interest area following the steps discussed in section "3. Manually Creating Interest Areas". In the property table of the interest area instance, set the "Left" value to 674, "Top" to 546, "Right" to 774, "Bottom" to 646, "Start Time relative to IP" to 0, and "End Time relative to IP" to 799. The subsequent interest area instances can be created using one of the methods described below.

Method 1: Manual Creation in Animation View

Set the trial view window to the animation view, as this view is useful for displaying, creating, and editing of instances of a dynamic interest area. You may play back the trial to the intended interest period time and manually draw each of the subsequent instances (please note that the top-left corner of the animation view displays both the current playback time relative to the currently selected interest period and to the start time of the trial). The coordinate, start time and end time can be adjusted through the property table of the instances.

Method 2: Copy and Paste

First copy the existing instance either from the event list in the Inspector Window or by clicking on it in the overlay / animation view and then paste it. If needed, for the newly created instance, edit the "left", "top", "right", "bottom", "Start Time relative to IP", and "End Time relative to IP" fields in the property table of the instance (the position of the interest area can also be adjusted by holding down the ALT key and then dragging it to the intended position). Repeat this process until all of the instances are created.

Method 3: Copy and Paste Current

First copy the existing interest area instance. Set the trial view window to the animation view and play back the trial to the intended interest period time (you may first drag the indicator on the slider at the top of the trial view window, and then use the left/right arrow key to get to the precise playback time). Now click the right mouse button and choose the "Paste Current" option. A new instance with proper time stamps will be created. You can adjust the position of the instance by holding down the ALT key and dragging the instance to its intended location, or by editing the "left", "top", "right",

“bottom” values in the property table of the instance. Repeat this process until all of the instances are created.

Method 4: Mouse Creation mode in animation view

First copy the existing interest area instance. Click on the right mouse button and choose the “Mouse Creation” option. Set the trial view window to the animation view and play back the trial to the intended interest period time (you may first drag the indicator on the slider at the top of the trial view window, and then use the left/right arrow key to get to the precise playback time). Now place the mouse cursor on top of the intended location and click the left mouse button – a new instance will be created following the mouse click. Repeat this process until all of the instances are created.

It may take some time for you to create the interest area instances for the very first recording file. Once the interest areas are saved into an .ias file, you can import that .ias file into the viewing session and use it as a template for subsequent similar trials across recording sessions.

5.8.25 Output Reports and Preferences

Version 2.0 of EyeLink Data Viewer supports report creation for both static and dynamic interest areas. However, dynamic interest areas introduce a number of complexities to data analysis. For example, you may run into scenarios where a single fixation lands in multiple interest areas. Some Data Viewer variables (e.g., the "CURRENT_FIX_INTEREST_AREA_INDEX" variable in the Fixation Report) require a fixation to be assigned to a single interest area. If a fixation lands in multiple dynamic interest areas, by default, this assignment will be based on the duration of fixation on respective interest areas - the interest area with the longest duration gets the fixation assignment. Occasionally, users may want to do the fixation assignment based on the index of the interest areas. This can be configured through the preference setting “Preferences -> Output/analysis -> Assign fixations to dynamic IA ordered by relative duration”. The default value is true, so to assign fixations by interest area index, uncheck this option to set it to false.

Another related issue when a fixation lands in multiple interest areas is the calculation of actual dwell time on each of the interest areas and fixation count. Data Viewer has an “Include Fixation in Single IA” option in “Preferences -> Output/analysis”. When selected, the fixation is assigned to one of the interest areas only (see the above paragraph on how the “winning” interest area is determined) - the interest area assigned gets the full length of the fixation when calculating dwell times. If this option is disabled (default setting), the fixation can be assigned to multiple interest areas, and effective duration (the total amount of time that fixation and instances of the same IA intersect) for each of the interest areas is used for the calculation of dwell times. This option will affect the calculation of some variables in the Fixation Report (e.g., CURRENT_FIX_INTEREST_AREAS, CURRENT_FIX_INTEREST_AREA_DWELL_TIME, etc.) and Interest Area Report (e.g., IA_DWELL_TIME, IA_FIXATION_COUNT, etc.). The following table illustrates

the effect of this option with four consecutive fixations. Columns “IA_1 Duration”, “IA_2 Duration”, “IA_3 Duration” are the duration of fixation onto each of three interest areas.

Fixation	Fixation Duration	IA_1 Duration	IA_2 Duration	IA_3 Duration	Winning IA
1	516	376	141	516	3
2	33	33	0	0	1
3	137	137	0	0	1
4	552	411	143	0	1

If the “Include Fixation in Single IA” option is checked, the first fixation is assigned to interest area 3, and the other three fixations are assigned to interest area 1. The IA_DWELL_TIME for interest area 1 will be 722 ms - the sum of full duration of the three fixations that are assigned to the interest area (33, 137, 552). The IA_DWELL_TIME value will be 0 ms for interest areas 2, and 516 ms for interest area 3. The IA_FIXATION_COUNT for each of the three interest areas will be 3, 0, and 1, respectively. If the “Include Fixation in Single IA” option is not checked, then the IA_DWELL_TIME calculation will be based on the effective duration (instead of the full fixation duration) on the target interest area. Thus, the IA_DWELL_TIME value for interest area 1 will be 957 ms ($=376 + 33 + 137 + 411$). The value for interest area 2 will be 286 ms ($=141 + 143$) and for interest area 3 it will be 516 ms. The IA_FIXATION_COUNT for each of the three interest areas will be 4, 2, and 1, respectively.

6 Data Analysis and Output

One of the ultimate goals of the EyeLink Data Viewer is to let you extract useful eye-movement information and export data to a tab-delimited text file that can be processed by most statistical packages. Before performing any data analysis, the first question you may ask yourself is what part of the recording data you need to look at? Do you need to analyze the eye movement data for the entire recording trial or just a small portion of it? For example, you conducted a reading experiment in which each trial displays a sentence followed by a comprehension question for the sentence just read. Do you need to look at the portion of eye movement data when the subject was answering the comprehension question? If not, you may create an interest period for the reading part (i.e., from the onset of the reading screen to the onset of the comprehension questions) and then select the intended interest period. In some experiments, you may create a reaction time period. The reaction time definition is a special variant of interest period filter in the sense that it reports the duration and end-event data in the Trial Report in addition to the data filtering capability.

Next, you will need to determine which report is appropriate for your analysis. EyeLink Data Viewer organizes the output reports based on different analysis emphasis. If you care about the summary statistics for a recording trial (e.g., how long is the trial recording, how many fixations were made in each of the trials?), the Trial Report is ideal for it. The Fixation Report, Saccade Report, and Sample Report provide detailed eye movement measures for each of the fixation, saccade events, or the individual eye samples. The Interest Area Report provides a columnar output for the eye-movement data related to each of the interest areas. The message report provides an output for each of the message events recorded in the data file with the corresponding fixation/saccade event. While most of the research questions can be answered by extracting and analyzing the data from one or several of those reports, it is possible that in some cases you will need to write some extra analysis code to further process the report outputs.

When performing analysis, you may organize each of the viewing sessions by loading one or multiple EDF files. The settings from one viewing session (a "template") can be exported to different viewing sessions. Thus the user may go through each of the output reports to configure the selection of variables in an intended order. Export the variable selection into a property file by clicking on the "Export Variable Selection" button. For all future viewing sessions, the variable selection in the reports can be automated by importing the previously saved property file (by clicking on the "Import Variable Selection" button). The export/import feature is similarly available to the interest period definitions and properties of the viewing session.

6.1 *Managing Trial Variables*

The Trial Variable Manager (Figure 6-1) and the Trial Variable Value Editor (Figure 6-2) allow the user to review the currently available variables, to assign a default value for the variables, to create or remove variables, and to change the variable values on a trial-by-trial basis.

6.1.1 Trial Variable Manager

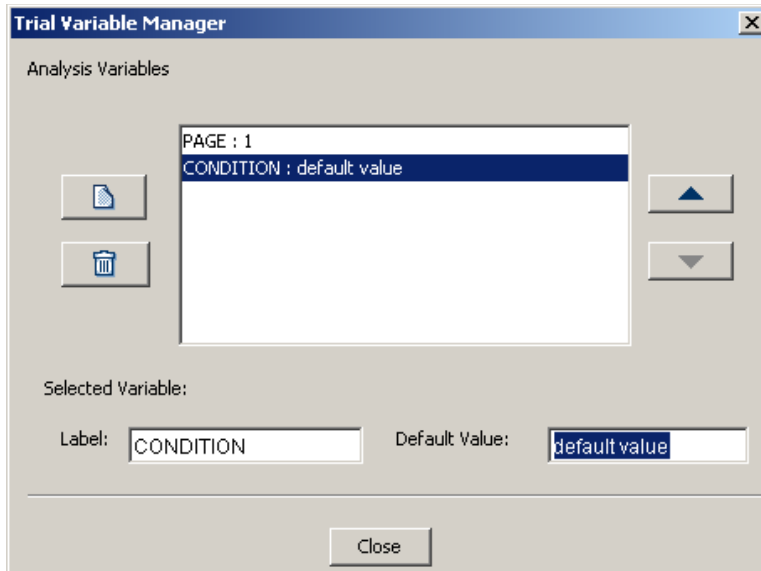






Figure 6-1. Trial Variable Manager


The Trial Variable Manager Dialog, activated by pressing **Analysis → Trial Variable Manager** from the menus, lists the filename (DATA_FILE) and other variables imported from the EDF file, if the TRIAL_VAR_LABELS command is used (PAGE in this case).

To assign a default value to one variable, click on the variable and type in the default value in the Default Value edit box (Be sure to press Enter to register the change you made). To create a new variable, click the  button on the left. A new variable (“label”) with a default value (“default value”) will appear. Select the variable and change its label and default value in the two edit boxes (DO NOT include quotes for text strings). To remove one variable, select it in the list and click the delete button .

Sometimes, the order in which the variables are listed is important. This is especially true when you want to match the variables created with the Trial Variable Manager and the values read from the command messages in EDF file (e.g., TRIAL_VAR_DATA, TRIALID; see the preference settings on data loading). To do that, simply select the variable and use the  or  button to move it to the desired position.

6.1.2 Trial Variable Value Editor

The variables and their values can be viewed and edited in the Trial Variable Value Editor dialog box.

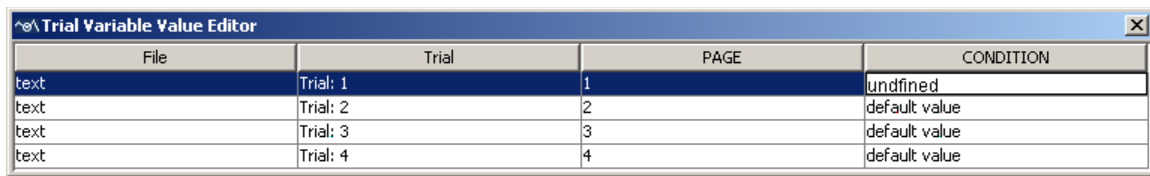

File	Trial	PAGE	CONDITION
text	Trial: 1	1	undefined
text	Trial: 2	2	default value
text	Trial: 3	3	default value
text	Trial: 4	4	default value

Figure 6-2. Trial Variable Value Editor Dialog

To activate the dialog box, from the menus, choose:

Analysis → Trial Variable Value Editor

The top-row of the dialog box lists the filename, variables imported from the EDF file, as well as variables created by the Trial Variable Manager. The following rows list the corresponding values for the variables in each individual trial. To change the value for a particular cell, double click on the current value, type in the new value (DO NOT include quotes for text strings), and press Enter to register the change. (Note, presumably, the user would load EDF files of the same experiment with the same condition labels, into one viewing session. Loading EDF files containing different condition labels from each other may result in some unexpected findings.)

Version 1.11.1 adds support for copying and pasting directly into the trial variable editor. If you need to add a new column into the editor, prepare an Excel sheet with the first row being the label of the column. Copy the selection from Excel and paste it onto top-left cell in the editor (the first row of the “File” column). If you already have the columns created but would like to update the content of the column, you can paste the content directly onto the first cell of the particular column without a row for the column label.

6.1.3 Trial Grouping

Trial condition variables are important for identifying one specific trial or grouping trials tested under the same experimental condition. Thus, instead of going through each of the individual trials, a trial grouping mechanism can be applied to manipulate those trials collectively. For example, practice trials can be removed from the viewing session in a batch, interest area templates can be applied to those trials using the same search display, and potentially, a “landscape” view can be created for the same display image with data collected from several subjects to identify the informative parts of the display.

By default, the Data Viewer group all of the trial by data file. To regroup trials within a viewing session, select the viewing session icon (✗), click the right mouse button, and choose the “Trial Grouping” option. In the following “Edit Trial Grouping” dialog, choose the variable with which the grouping should be performed and press the

“Regroup” button (see the left panel of Figure 6-3). Multiple grouping variables can be selected or de-selected by using the control key; if no grouping variable is chosen, grouping will be performed based on the data file each trial belongs to. Once grouping is performed, the trials will be listed by the conditions of the grouping variables selected (see the right panel of Figure 6-3). The property panel of each group lists the label (values of the grouping variables), total number of trials, and configurable interest area information for the group. Thus, the user can potentially apply an interest area template to all trials in the group (see section 5.8.7 Using Interest Area Templates).

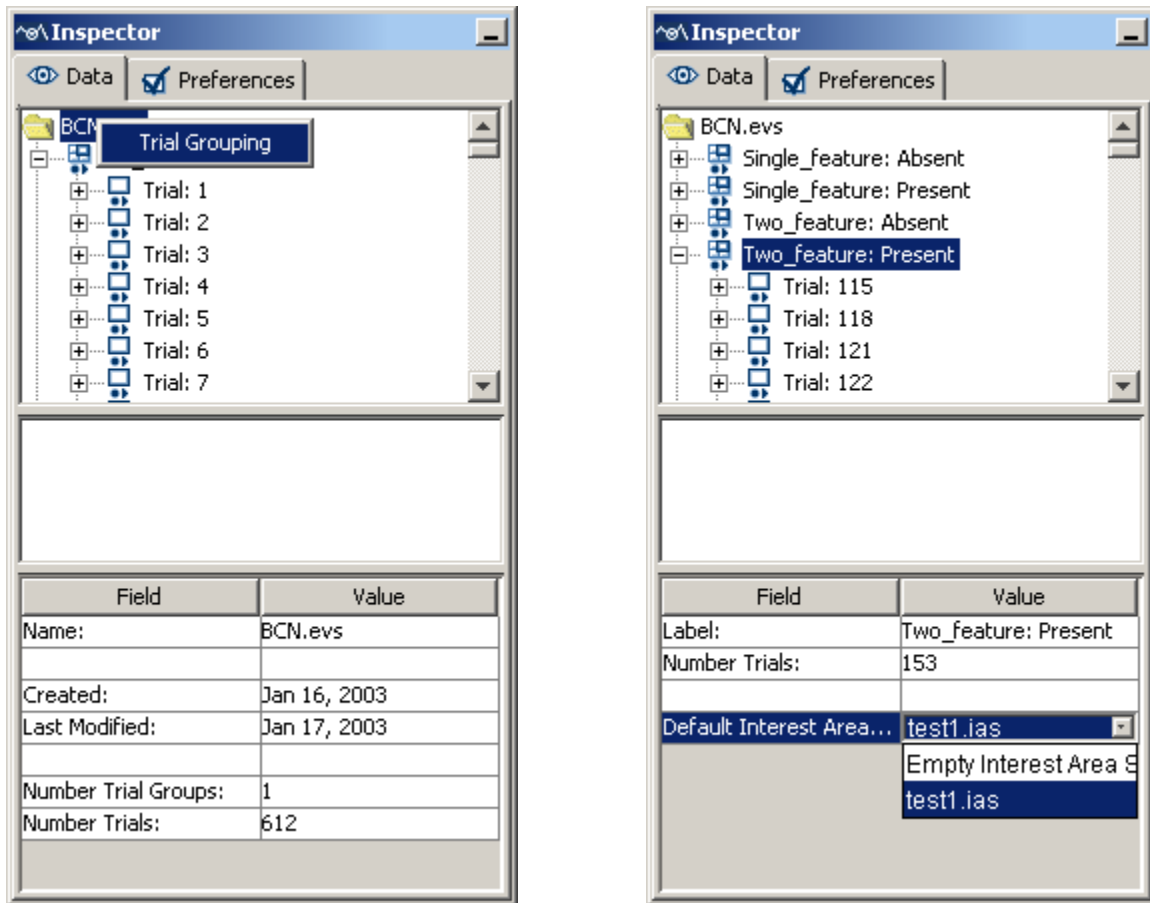


Figure 6-3. Applying Trial Grouping by a Trial Condition Variable and Setting Interest Area Template for a Group of Trials.

Version 1.11.1 adds a “Report Hidden” option for the groups so that users can block outputting particular groups of trials by enabling this option.

6.2 Fixation Map

EyeLink Data Viewer allows users to create a "landscape" view to identify informative parts of a display based on recording trials from a single participant or multiple participants with the same background image. To create a fixation map (or heat map), a

2D Gaussian is applied to each of the fixation - the Gaussian center is the fixation location, the width of the Gaussian is influenced by an adjustable sigma value in degrees of visual angle (the larger the sigma value, the larger the area that will be influenced by the fixation), and the height of the Gaussian can be weighted either equally for all fixations or by the duration of individual fixations. This 2D Gaussian is added to an internal map by adding weight to that area of the map. After the above process is applied for all fixations, the internal map is normalized so it can be applied to the color or brightness scale used to create an aggregate view (see Figure 6-4).



Figure 6-4. Example fixation map

To create a fixation map for a single trial, select the trial node (🖥️) in the top-panel of the inspector window, click the right mouse button and select "Create Fixation Map" option (see Figure 6-5). To create a fixation map for a group of trials, select the group node (📁). Click the right mouse button and select "Create Fixation Map" option (if no common image has been applied to the trials yet, select "Select Trial Background Image" option before creating a fixation map). In the following "Save Fixation Map ..." dialog box, enter the intended file name with a proper file extension .BMP, .JPG, or .TIFF (a .bmp file will be created by default if no file extension is specified). Properties for a fixation map can be configured through the Output/Analysis preferences settings.

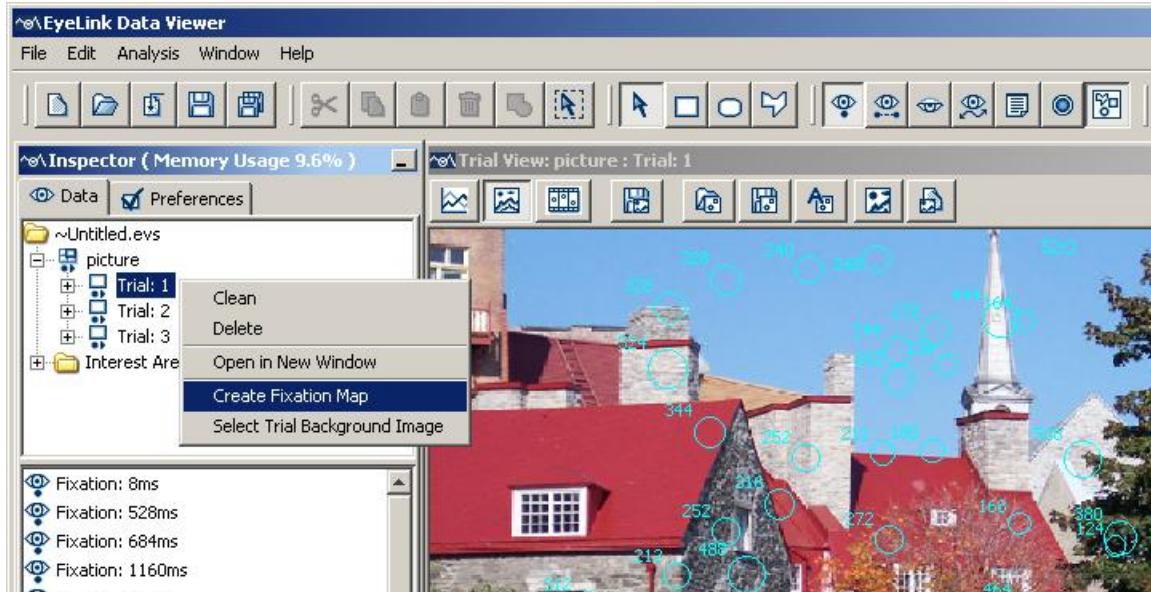
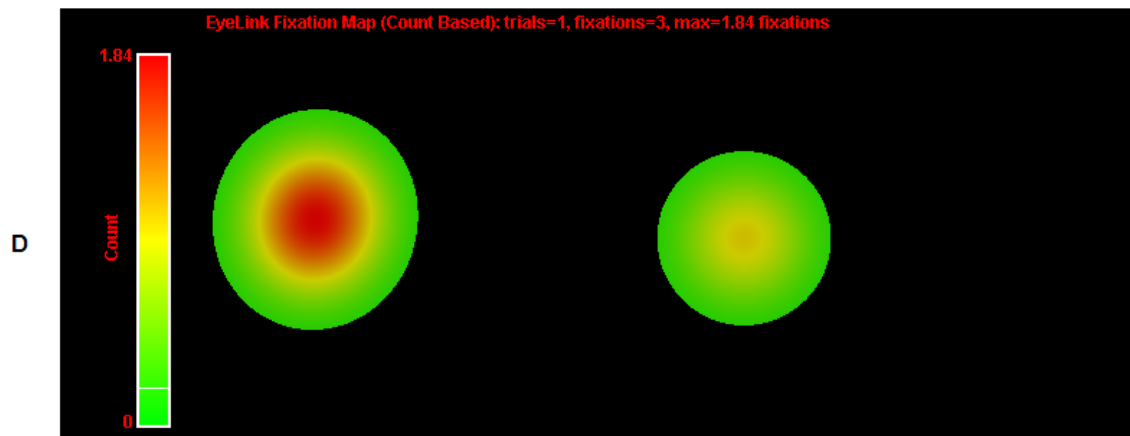
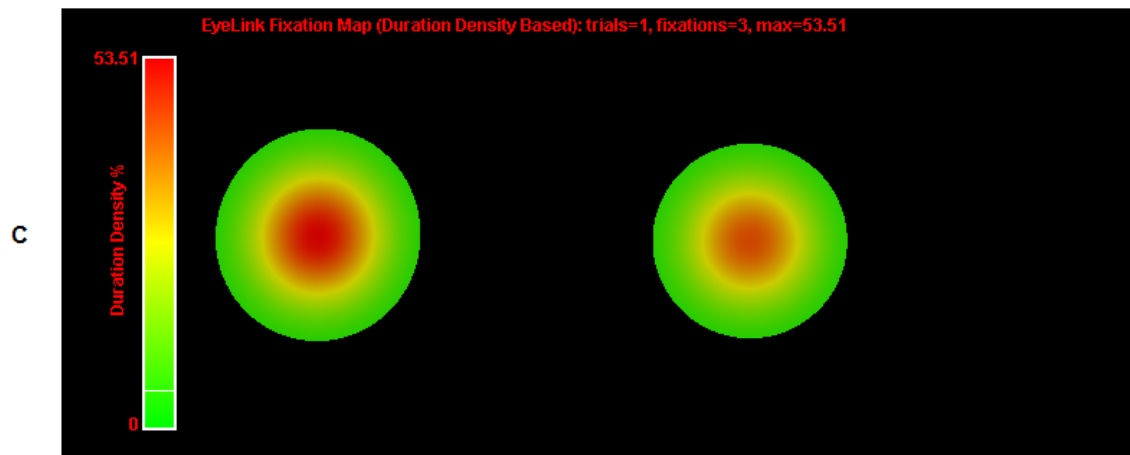
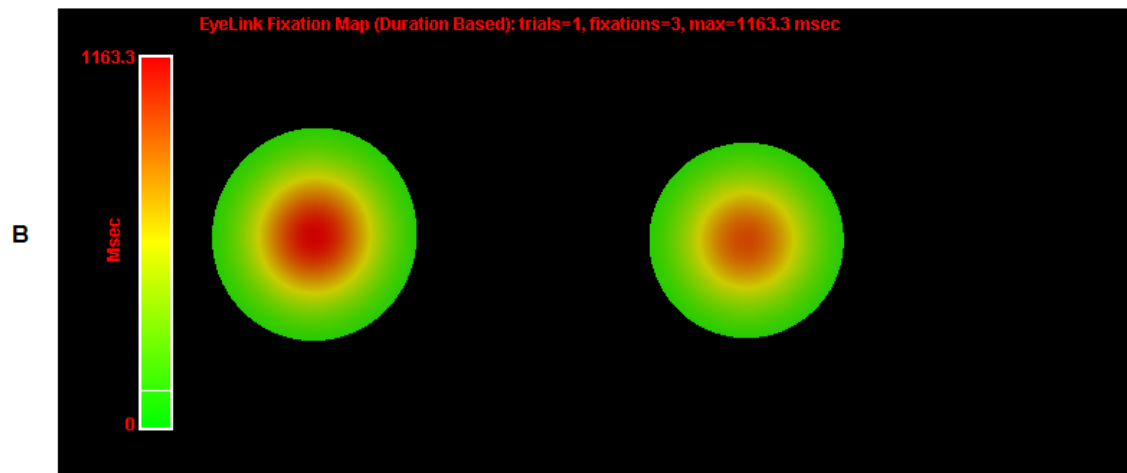


Figure 6-5. Creating a Fixation Map

EyeLink Data Viewer supports creating a fixation map based on the duration of fixations, count of fixations, percent of fixations, and percent of dwell time across the display. This is configurable through the "Type (1 = Duration, 2= Count, 3 = Count Density, 4 = Duration Density)" option of the Output/Analysis preferences settings. If a duration or duration-density type is selected, the height of the Gaussian is weighted by the duration of the individual fixations; for a count-based or a count density fixation map, the height of the Gaussian is the same for every fixation. The map activity is averaged based on trial count pooled across subjects (i.e., a map based on 1 trial that has 1 fixation will have a same fixation output as a map based on 10 trials, each trial with 1 fixation at same location). The title of the fixation map indicates the type of fixation map created, the total number of trials and fixations contributing to the output, and the maximum per-trial-average value for the type of map chosen. The legend of the map indicates how long or how often subjects have looked across the map. By default, the frequently fixated areas are painted in red whereas the less-fixated areas are painted in green. The drawing colors and the transparency of the drawing can be configured through the preferences. A gray-scale version is also available if the "Enable Heat Map Mode" option is unchecked.

The following is a sample output of the four different fixation maps for the same data trial: A - original fixations in the spatial overlay view; B - the duration-based fixation map; C - the density distribution of the fixation durations (i. e., proportion of trial dwell time on each area of the display); D - the count-based fixation map; and E - the density distribution of the fixation counts (i.e., how frequently each area of display area has been visited).





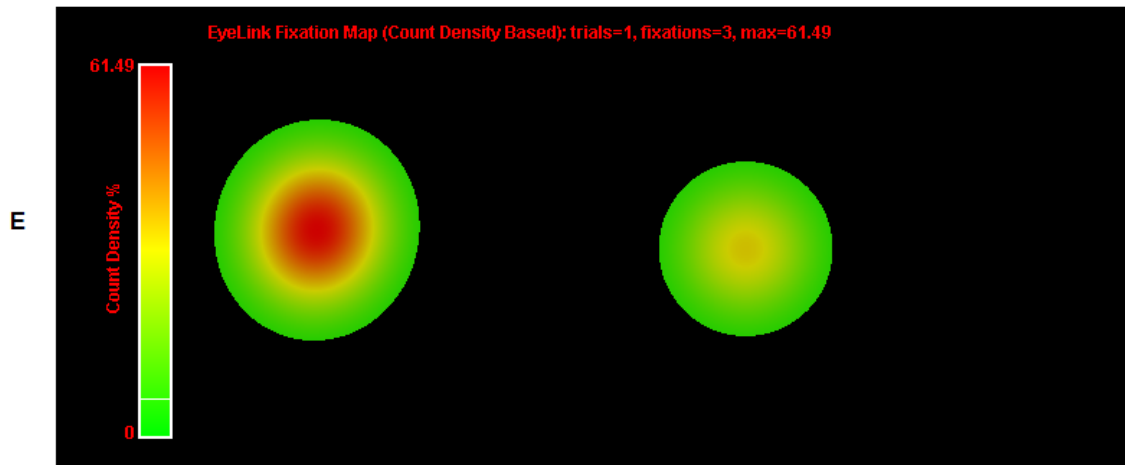


Figure 6-6. Types of Fixation Map

A single fixation map is useful for the purpose of identifying informative parts of a display; the legend displays colors that correspond to the peak and lowest activation values. However, fixation activities across two different fixation maps cannot be directly compared if the scales of the two maps are very different. To make such a comparison, the user may set an intended peak activity value in the "Fixed Maximum Value" field of the preference settings. By default (with a value of -1 in the "Fixed Maximum Value" field), the peak value displayed in the legend will be the actual maximum per-trial-average value for the map. If a user-defined peak value is set, the drawing of fixation map will be adjusted according to this new scale. A black line is drawn on the heatmap scale to indicate the actual maximum activation of the heatmap (if that value is less than the fixed maximum). A white line is drawn on the heatmap scale to indicate the "Low Activity % Cut-off" when the setting is greater than 0 and the "Fixed Minimum Value" set in the preference. This low activity cutoff feature ignores areas on the map with an activation value lower than 10% (default) of the peak value. If the low activity cutoff value is set to 0, the whole display is painted.

Panel A of the following figure shows the original fixations in a trial. Panel B shows the fixation-count based map using the default "Fixed Maximum Value" (-1) and "Low Activity % Cut-off" (10% or 0.2 fixation - the white line position). Panel C shows the same trial but with a different "Fixed Maximum Value" (4) - the actual peak value in the map is indicated by a black line. Panel D illustrates the output using a "Fixed Maximum Value" of 2.0 (which is the same as the actual peak value) and a "Fix Minimum Value" of 0.5 (which is higher than the default 10% cutoff value, i.e., 0.2 fixation) with areas below the fixed minimum value clipped.



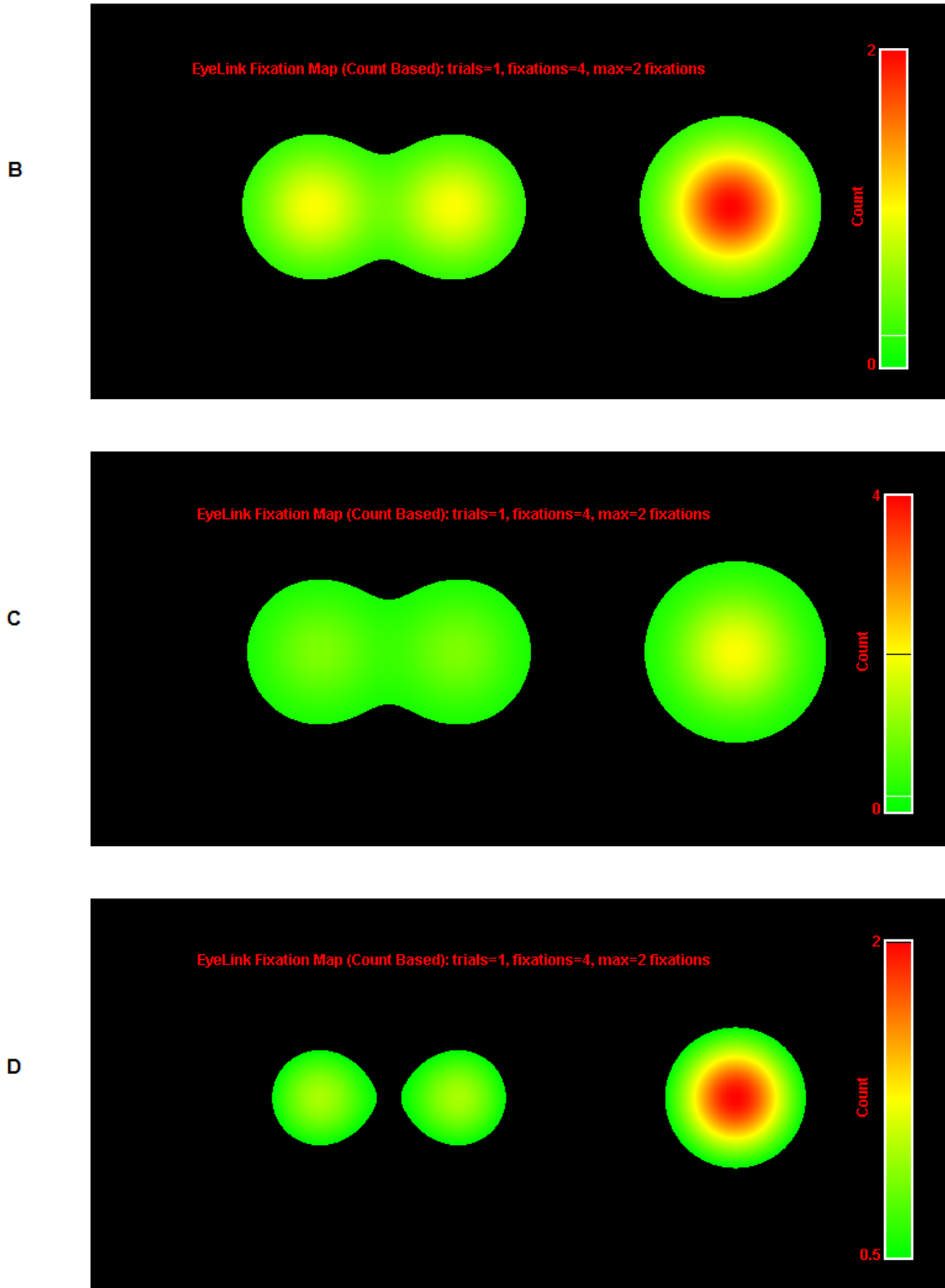


Figure 6-7. Adjusting the Scale of the Fixation Map

6.3 Managing Reaction Time Definitions

Another important application of the trial condition variables is for managing reaction-time definitions for all the trials in the same experimental condition. A typical trial recording could be ended in several ways, such as a button response, a saccade event, a fixation event, or a special message. To obtain an accurate reaction time (RT) measure, the EyeLink Data Viewer relies on a set of reaction time definitions to parse the start and end events for RT calculation. The reaction time manager allows the user to configure the way in which the trial starts and ends for each condition of the experimental design. Once a set of RT definitions has been created and initialized, the RT start and end event information is available in both the trial view window and the inspector window.

6.3.1 Reaction Time Manager

The Reaction Time Manager is used to create, edit, or delete reaction time definitions for different experimental conditions. The user can set parameters of a trial end event and specify condition values that must be matched for the RT definition to be applied to a given trial.

To configure reaction time definition,

- 1) From the menus, choose:
Analysis → Reaction Time Manager

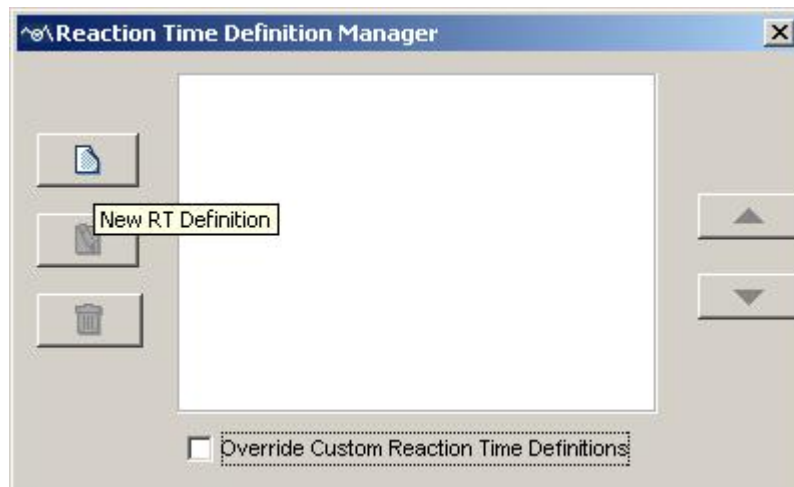






Figure 6-8. Reaction Time Definition Manager

In the reaction time manager dialog (see Figure 6-8), the user can create a new reaction time (RT) definition by clicking the new  button on the left, edit an existing RT definition by clicking on the  button, and remove an RT definition by clicking on the  button.

Users who have already had custom reaction time messages defined in the EDF file need to decide whether the customer RT messages or the RT definitions

created here should take precedence when both of them are applicable for a particular trial. If "Override Custom Reaction time definitions" box is checked, the new RT definitions created through the Reaction Time Manager will take precedence; otherwise, the custom RT messages written in the EDF file will be used.

- 2) Click on the  button. In the Reaction Time Definition Editor dialog (General Tab; see Figure 6-9)
 - a. Enter a label for the current RT definition in the Label edit box;
 - b. Select the type of RT end event from the Type combo box;
 - c. In the Trial Condition Qualifiers group box, select the trial condition variables and enter the intended values for the current RT definition (DO NOT include quotes for text strings).

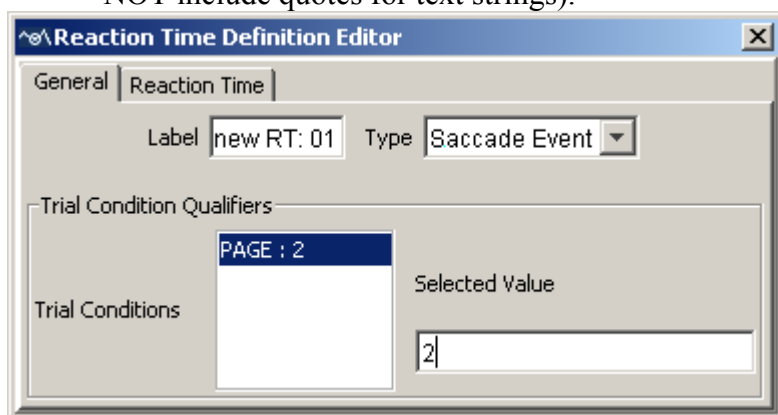


Figure 6-9. Reaction Time Definition Editor (General Tab)

- 3) Select the Reaction Time tab (see Figure 6-10).

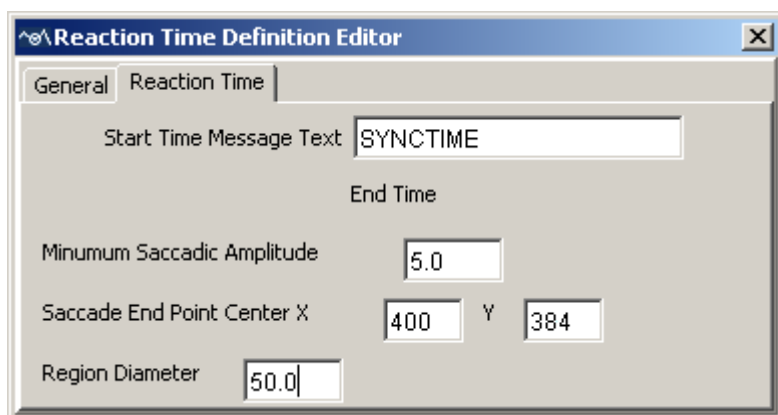


Figure 6-10. Reaction Time Definition Editor (Reaction Time Tab)

The layout of the Reaction Time tab depends on the type of end event you selected in the General tab. One common feature though, is the configuration of

the start time message. In the Start Time Message Text edit box, type in the intended message string (the default is "SYNCTIME"). The "Use Substring Match for Messages" option, if checked (default), searches for any string that contains the message text specified here. If this option is unchecked, the message string should be exactly like the one specified. For example, "SYNCTIME_SINGAL" will not be accepted if the specified text is "SYCTIME" and the "Use Substring Match for Messages" option is unchecked.

Button Event:

This will treat a particular button press/release as the reaction time end event. Select the intended button ID from the Button Number combo box and choose the button state (button press or release).

Message Event:

This will treat the first occurrence of a matching message as the end event for reaction time calculation. In the "Message Text" edit box, type in the message string (DO NOT include quotes for text strings).

Fixation Event:

This will end the trial by the appearance of a specific fixation. In the edit boxes, fill in the minimum fixation duration (in millisecond) required, as well as the *xy* coordinates of the center of the fixation region and the diameter of the region.

Saccade Event:

This will end the trial by the occurrence of a specific saccade. Fill in the minimum saccadic amplitude (in degrees of visual angle) required, as well as the *xy* coordinates of the region center and the diameter of the saccade region.

- 4) Once all the desired RT definitions have been set, arrange the order of the definitions by using the ▲ and ▼ buttons in the Reaction Time Definition Manager dialog. This operation is important only when more than one RT definition can be met for a single trial. Always put specific definitions before general ones.
- 5) Close the Reaction Time Manager dialog. This will apply the reaction definitions to all trials. If there are lots of trials in a viewing session, this may take some time to finish.
- 6) If the user thinks that the current RT definitions are not appropriate, she/he can remove them by clicking on the "Delete" button on the Reaction Time Definition Manager dialog box.

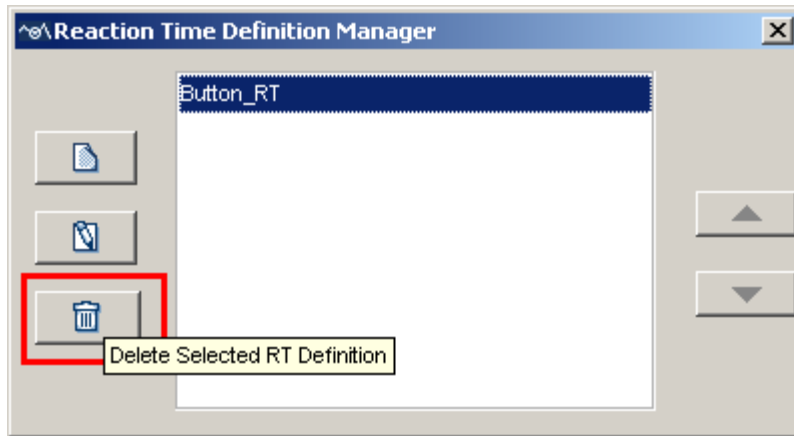



Figure 6-11. Deleting Reaction Time Definitions

6.3.2 Viewing Reaction Time Information

If the reaction time events are initialized, such information is available for viewing in trial view window (see Figure 6-12). In the spatial overlay view, a red  symbol is displayed near the end event. In the temporal graph view, a green line is drawn at the time when the start time message is presented and a red line is drawn at the time when the end event occurs (to see these graphics, make sure that the Display RT Graphics field in the General Trial View Preferences is checked; see section 8.5).

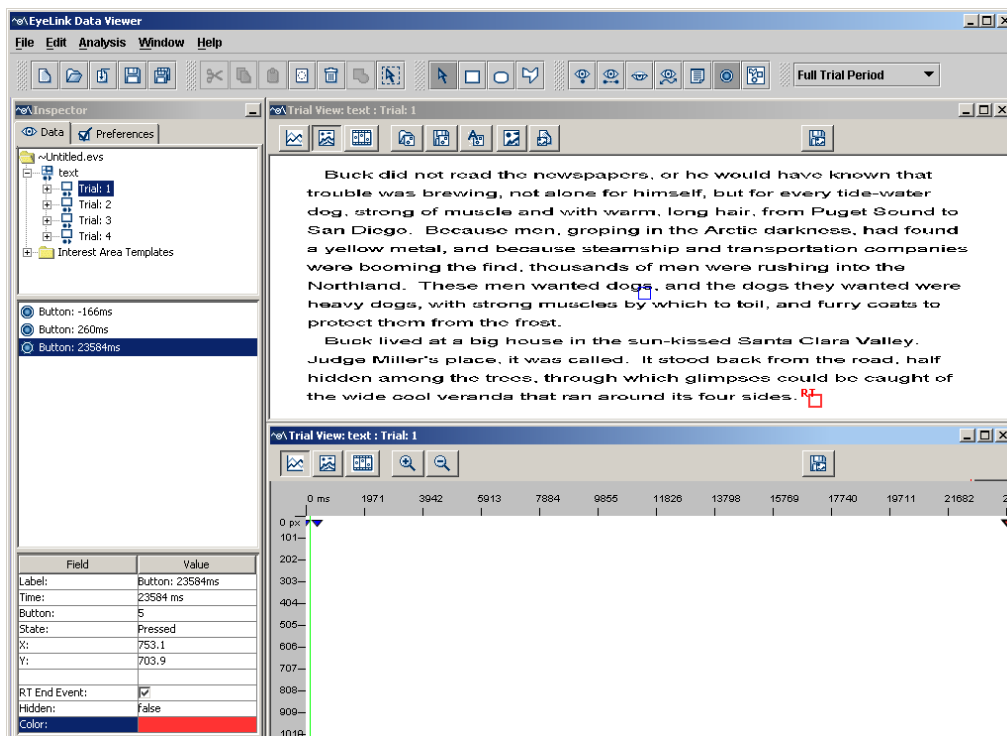


Figure 6-12. Reaction Time Graphics in Both the Spatial Overlay View and the Temporal Graph View of the Trial View Window






In the event list of the Inspector window, select the end event that meets the reaction time definition for the selected trial. The RT End Event field of that event is automatically checked. This field will be unchecked and reaction time calculation will be updated if the user decides to choose another event as the RT end event by checking the same field of the newly selected event.


6.4 Interest Periods

Besides the above-mentioned reaction-time definitions, the user may also selectively view and perform analysis on data within a specific period of time within a trial (Interest Period). For example, the user may want to examine the eye data between two important markup messages; for a trial with multi-page manipulation, analysis may be performed separately for the period of each page presentation. The EyeLink Data viewer allows the user to create interest periods based on messages, button events, or duration. To do that, when programming an experiment, the user should mark up possible interest periods with messages or other important events such as button presses. If different images were shown within different interest period, the user should also make sure that messages used to load images, text, or draw simple graphics be recorded with the scope of the intended interest period.

6.4.1 Creating Interest Periods

To create an interest period,

- 1) Click the "Interest Period List" combo box at the far-right hand of the application menu bar and select the "Edit ...".
- 2) Similar to the reaction time manager dialog, the user can create a new interest period by clicking the New  button on the left, edit an existing interest period by clicking on the Edit  button, and remove an interest period by clicking on the Delete  button. Data Viewer also supports exporting and importing of interest periods across viewing sessions. For example, the user may create intended interest periods for one session. Export the interest periods into a property file by clicking on the "Export the Interest Period" button . For all future viewing sessions, creation of the interest periods can be automated by importing the previously saved property file (by clicking on the "Import the interest period" button ).

Click on the new  button. In the Interest Period Editor dialog (General Tab, see Figure 6-13), enter a label for the current interest period in the label edit box and select the type of interest period end event (Button, Message, or Duration-based) from the Type combo box.

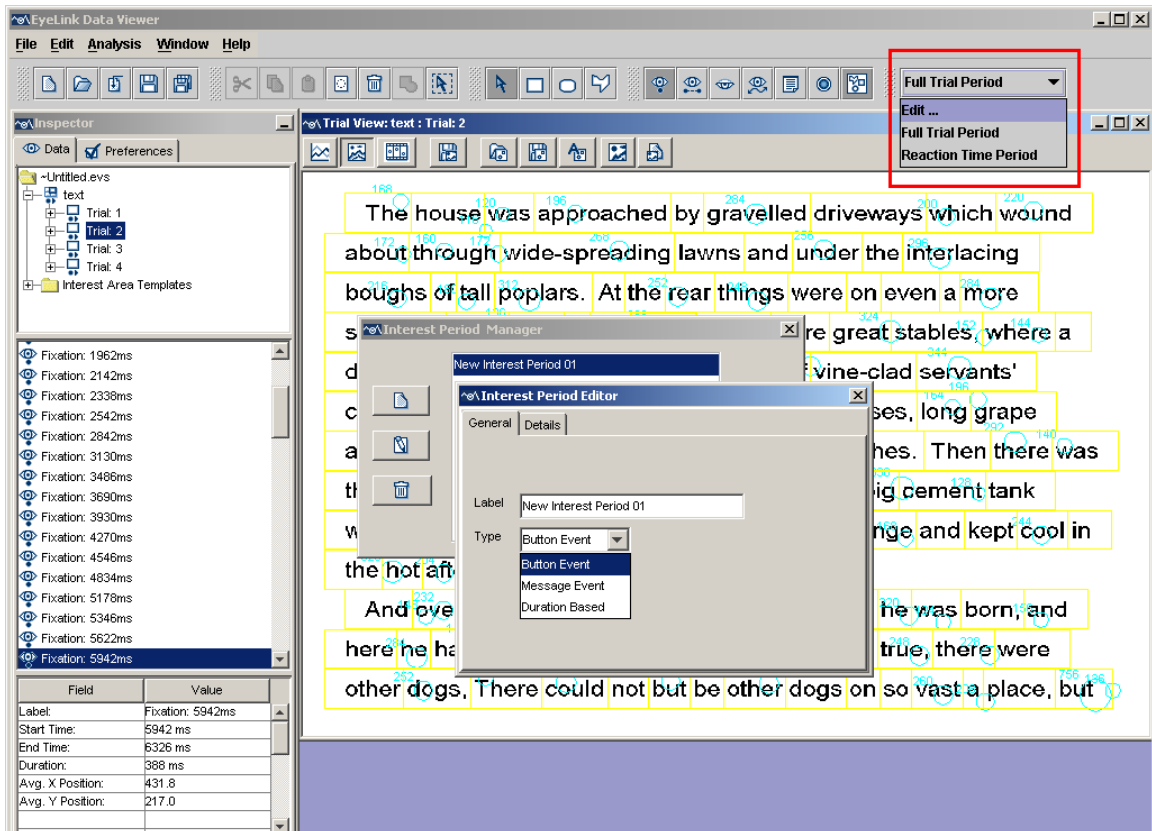


Figure 6-13. Interest Period Editor (General Tab)

3) Select the Details Tab (see Figure 6-14).

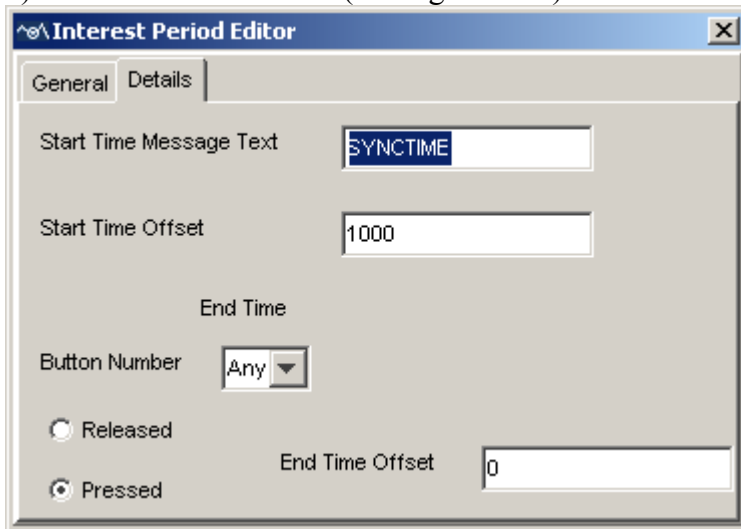


Figure 6-14. Interest Period Editor (Details Tab).

Similar to the reaction time definition, the layout of the Details tab depends on the type of the end event the user selected in the General Tab. A "Use Substring Match for Messages" option is available for the "Start Time Message Text" and "End Time

Message Text". If checked (default), Data Viewer searches for any string that contains the message text specified here. If this option is unchecked, the message string should be exactly like the one specified. For example, "SYNCTIME_SINGAL" will not be accepted if the specified text is "SYCTIME" and the "Use Substring Match for Messages" option is unchecked.

Button Event:

This will treat a particular button press/release as the interest period end event. In the Start Time Message Text edit box, enter the message string (case-sensitive; DO NOT include quotes for text strings); if an empty string is used, this will be the start of the trial recording. Select the intended button ID from the Button Number combo box and choose the desired button state (button pressed or released). The user can also specify a time offset (0 or a positive number in milliseconds) for the start time and end time so that the interest period does not start or end right on the time of message and/or button event.

Message Event:

This will treat the first occurrence of a matching message as the end event for the interest period. Enter the desired message strings for both the Start Time and End Time Message Text edit boxes (case sensitive, with no quotes for text strings); a valid message must be supplied for both start and end time message text fields. The user can also set a time offset (0 or a positive number, in milliseconds) for the start time and end time. Please note that when using this option, the interest period end message must be different from the start message; otherwise, the user should use a duration-based interest period instead (see below).

Duration Based:

The interest period can also be based on a duration following the detection of the start message (e.g., 10 seconds following the "SYNCTIME" message). The user needs to specify a message string (case-sensitive; DO NOT include quotes for text strings) in the Start Time Message Text edit box; if an empty string is used, this will be the start of the trial recording. The user can also specify an optional offset time value so that the interest period does not start right on the message time. In the "Duration" edit box, enter the desired duration (a positive number in milliseconds) of the interest period.

4) Close the Interest Period Editor dialog. This will initialize the interest period for all trials.

6.4.2 Using Interest Period

The newly-created interest periods are added to the interest period list, along with the default interest periods – "Full Trial Period" and "Reaction Time Period". The user can navigate between different interest periods by selecting the desired interest period from the list. For each interest period, only those events falling within that period are shown in the spatial overlay, temporal graph views, and in the event list of the inspector window.

In addition, for all of the following output analyses (trial, fixation, saccade, interest and recording sequence outputs), only those events and data within the selected interest period (custom-defined interest period, or the default – “Full Trial Period” or “Reaction Time Period”) are exported.

6.5 Trial Report

Trial report summarizes the performance of individual trials. Before trying to obtain a trial report, make sure that you have completed the reaction time definition and initialized the reaction time events.

6.5.1 Obtaining a Trial Report

To obtain a trial report,

- 1) From the application menus, choose:
Analysis -> Report -> Trial Report

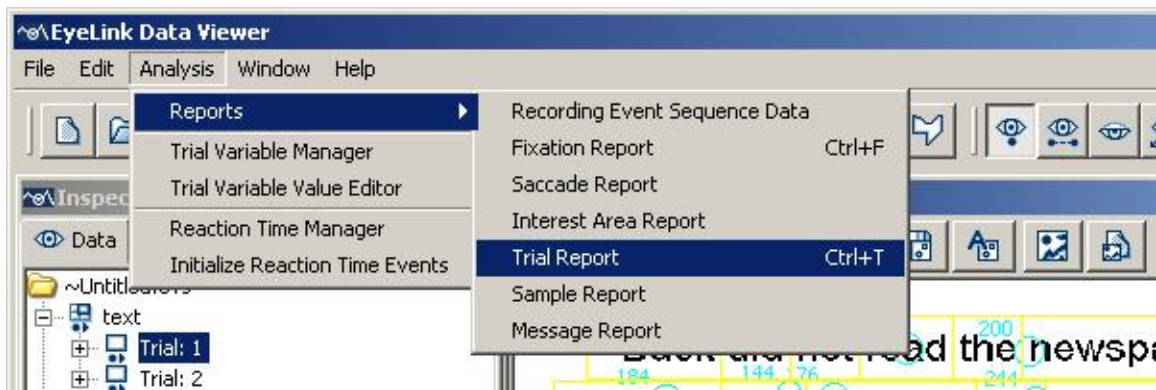


Figure 6-15. Creating a Trial Report.

- 2) In the following output report dialog, select the variables in the list of available output variables (left panel) and press ">>" to enter the variables into the list of selected output variables.

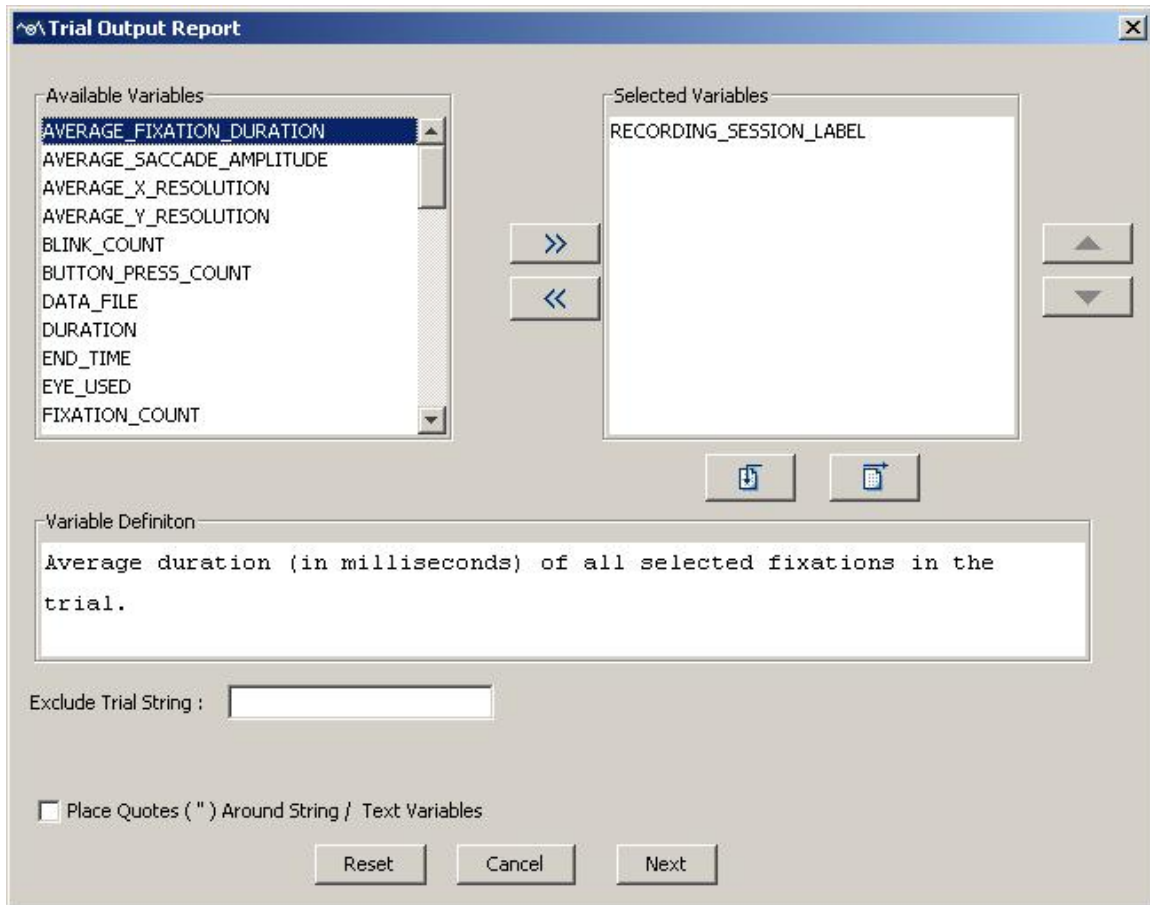




Figure 6-16. Variables used in the Trial Report.

- To remove one variable from the output list, simply select that variable and press <<.
- To change the order in which the variables are listed, simply select one variable and use the ▲ or ▼ button to move it to the desired position. Repeat this step until all variables are in the right position.
- Data Viewer supports exporting and importing of variable selection across viewing sessions. For example, the user may go through each of the output reports to configure the selection of variables in an intended order. Export the variable selection into a property file by clicking on the "Export Variable Selection" button . For all future viewing sessions, the variable selection in the reports can be automated by importing the previously saved property file (by clicking on the "Import Variable Selection" button .

3) Once all the desired variables have been entered into the output list, press the Next button.

4) In the Export dialog box, browse to the directory in which you want to save the output report.

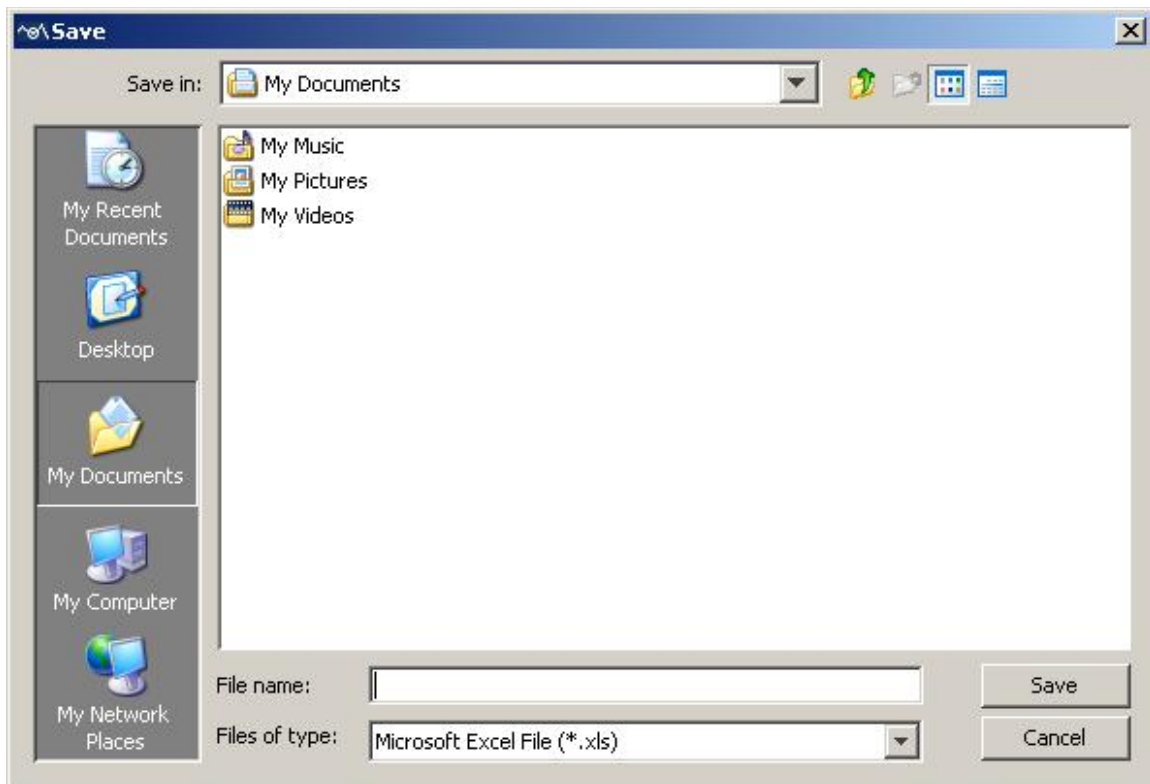


Figure 6-17. Saving the Output Report.

5) In the File name text box, type in the name for the document and choose the appropriate file extension. Note, if the output report file extension set to "All Files" and the file name does not contain an extension, the file name is appended with a ".txt" by default.

6) Press the "Save" button.

6.5.2 Variables in the Trial Report

The following variables are listed in the trial report:

Field	Contents
RECORDING_SESSION_LABEL	Label of the data file
* AVERAGE_BLINK_DURATION	Average duration (in milliseconds) of all selected blinks in the trial.
* AVERAGE_FIXATION_DURATION	Average duration (in milliseconds) of all selected fixations in the trial
* AVERAGE_SACCADE_AMPLITUDE	Average size (in degrees of visual angle) of all selected saccades in the trial
AVERAGE_X/Y_RESOLUTION	Average horizontal/vertical angular resolution

	(in screen pixels per degree) for the trial
* BLINK_COUNT	Total number of blinks in the trial
* BUTTON_PRESS_COUNT	Total number of button presses in the trial
DATA_FILE	File name of the data
DURATION	Duration of the trial recording between the “START” message and the “END” message
END_TIME	Timestamp when the trial recording ends (in milliseconds since EyeLink tracker was activated)
EYE_USED	Which eye's data (LEFT or RIGHT) is used to create this report?
* FIXATION_COUNT	Total number of fixations in the trial
FIXATION_DURATION_MAX	The longest fixation duration (in milliseconds) in the trial.
FIXATION_DURATION_MIN	The shortest fixation duration (in milliseconds) in the trial.
FIXATION_DURATION_MAX_TIME	Start time (in absolute EDF file time) of the fixation with the longest fixation duration in the trial.
FIXATION_DURATION_MIN_TIME	Start time (in absolute EDF file time) of the fixation with the shortest fixation duration in the trial.
IA_COUNT	Reports total number of unique interest areas in the trial.
INDEX	Sequential order of the trial in the recording
INTEREST_AREA_SET	Setting of the interest area (“Custom Interest Area Set” if interest areas are loaded/created and “Empty Interest Area Set” if not)
* IP_DURATION	Duration of the interest period in milliseconds.
* IP_END_TIME	End time (in milliseconds since EyeLink tracker was activated) of the interest period.
* IP_START_TIME	Start time (in milliseconds since EyeLink tracker was activated) of the interest period.
* MESSAGE_COUNT	Total number of messages in the trial
* PUPIL_SIZE_MAX	Largest pupil size in arbitrary units (Typical pupil area is 800 to 2000 units while pupil diameter is in the range of 1800-3000 units) in the trial recording. Note that PUPIL_SIZE_MAX, PUPIL_SIZE_MEAN, and PUPIL_SIZE_MIN fields will output a missing value if samples are not loaded into a

	viewing session (see Data Loading Preferences).
*PUPIL_SIZE_MAX_TIME	EDF Time of the sample with maximum pupil size. Note that this field will output a missing value if samples are not loaded into a viewing session (see Data Loading Preferences).
*PUPIL_SIZE_MAX_X	X position at PUPIL_SIZE_MAX_TIME.
*PUPIL_SIZE_MAX_Y	Y position at PUPIL_SIZE_MAX_TIME.
*PUPIL_SIZE_MEAN	Average pupil size in arbitrary units in the recording trial.
*PUPIL_SIZE_MIN	Smallest pupil size in arbitrary units in the recording trial.
*PUPIL_SIZE_MIN_TIME	EDF Time of the sample with minimum pupil size.
*PUPIL_SIZE_MIN_X	X position at PUPIL_SIZE_MIN_TIME. Note that this field will output a missing value if samples are not loaded into a viewing session (see Data Loading Preferences) or the eye is in a blink.
*PUPIL_SIZE_MIN_Y	Y position at PUPIL_SIZE_MIN_TIME. Note that this field will output a missing value if samples are not loaded into a viewing session (see Data Loading Preferences) or the eye is in a blink.
¶ REACTION_TIME	Configurable response time determined by the Reaction Time Definition associated with the trial. This variable returns a '.' if the trial does not have an RT end event.
¶ RT_DEFINITION_LABEL	Label of the RT definition (“.” if undefined). “CUSTOM_RT” if the RT definition is a custom one.
¶ RT_EVENT_BUTTON_ID	If the RT end event is the press/release of a button, records the button ID; otherwise, records “.”
¶ RT_EVENT_END_TIME	End time (in milliseconds since EyeLink tracker was activated) of the end event. If end event is not an eye event (fixation or saccade event), this variable is equal to RT_EVENT_START_TIME.
¶ RT_EVENT_INDEX	Sequential order of the end event; -1 when NO_MATCH. Index is relative to a given event type.

¶ RT_EVENT_MESSAGE_TEXT	Records the message text of the end event. If the end event is not a message event, records “.”
¶ RT_EVENT_POSITION_X/Y	Records the X/Y coordinates of the fixation or saccade end event and sets “.” for a button or message end event.
¶ RT_EVENT_START_TIME	Start time (in milliseconds since EyeLink tracker was activated) of the end event.
¶ RT_EVENT_TYPE	The way in which the trial is ended (BUTTON, FIXATION, SACCADE, MESSAGE, or NO_MATCH)
¶ RT_START_TIME	Start time (in milliseconds since EyeLink tracker was activated) of RT definition for a trial.
* RUN_COUNT	Total runs of fixations in the trial (two consecutive fixations in the same interest area belong to the same run).
* SACCADE_COUNT	Total number of saccades in the trial
* SAMPLE_COUNT	Total number of samples in the trial
START_TIME	Timestamp when the trial recording starts (in milliseconds since EyeLink tracker was activated)
VISITED_INTEREST_AREA_COUNT	Reports total number of unique interest areas visited in the trial (i.e., RUN_COUNT without replacement).
TRIAL_LABEL	Label of the trial

Besides these default variables, additional trial variables will be listed if they have been defined in the EDF file or created during the viewing session (see section 6.1.1 on trial variable manager).

6.5.3 Using Trial Report

Please note that in the above table, outputs for all of the variables marked with an “*” will change as different interest period filters are used. For example, the output for a trial filtered by a reaction time period may be different from a full-trial output, as the former only include a subset of the full-trial data in the report. Therefore, please make sure that you have the right setting for the interest period filter when creating you own reports (see figure below; see also 6.4.2). In those trials with no reaction time defined, missing values (“.”) will be recorded for those variables marked with a “¶” sign. Please also note that event counts only factor in visible events. If an event is hidden and not visible, it will not contribute to event counts. If hidden events are visible, then they will apply to event counts.

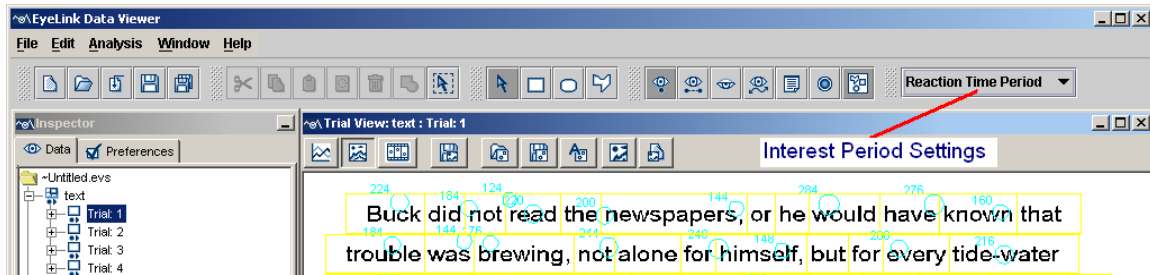


Figure 6-18. Using Interest Period for Event Filtering

Three time periods have been used in a trial output report: 1) recording period, 2) reaction-time period, and 3) interest period. The recording period (START_TIME, END_TIME, and DURATION) of a trial starts from the “START” message in the EDF file and ends on the “END” message. This is fixed for a recording trial, regardless how the user set the interest period filters.

If a valid reaction time definition is found for one trial, the RT_START_TIME is the time of the RT start message (e.g., time of “SYNCTIME” message), RT_EVENT_END_TIME, and RT_EVENT_START_TIME are the start and end time of the RT event (fixation, saccade, button, or message). Reaction time (REACTION_TIME) is calculated as the difference between RT_START_TIME and RT_EVENT_START_TIME. Please note that the RT_EVENT_END_TIME and RT_EVENT_START_TIME will be the same when a button or message is used as the RT end event but will be different when a fixation or saccade is defined as the end event. If the user does not create a reaction time definition or no valid reaction time definition is found for one trial, REACTION_TIME, RT_EVENT_END_TIME, and RT_EVENT_START_TIME, and other variables related to reaction-time definition will have missing values (“.”) whereas RT_START_TIME will be the same as the trial recording start time (START_TIME).

The interest period variables will have different values depending on the settings used in the interest period filter (see Figure 6-16). If the interest period filter is set to “Full Trial Period”, the start (IP_START_TIME), end (IP_END_TIME), and duration (IP_DURATION) of the interest period will be same as those for the recording period (i.e., START_TIME, END_TIME, and DURATION). If the interest period filter is set to “Reaction Time Period”, the start, end, and duration of the interest period will be same as those for the reaction time period. If the user has created her/his own interest period (see section 6.4.1) and that interest period is used to filter events, the output report will only do summary statistics (variables marked with “*” in the above table) for those events falling within that period.

6.5.4 Configuring the Trial Report

The Data Viewer also allows the user to configure the trial report (similarly the fixation report, saccade report, and the interest area report). The user can exclude some trials by entering a string in the “Exclude Trial String” edit box. This string should be one message or part of a message within the scope of the trial recording (see section 7.1 on defining the start and end of a trial) that uniquely defines that subset of trials to be excluded. In addition, to enable correct loading of string/text variables in some statistics software, the Viewer also has an option of adding a pair of quotation marks for such variables.

6.6 Fixation Report

Fixation report creates a columnar output of fixation events. Each row of the report represents a fixation event. Depending on the variables selected for output, each row can contain information on the previous fixation and/or saccade, as well as the next fixation and/or saccade.

6.6.1 Obtaining a Fixation Report

To obtain a fixation report ,

- 1) From the application menu, choose:
Analysis -> Report -> Fixation Report

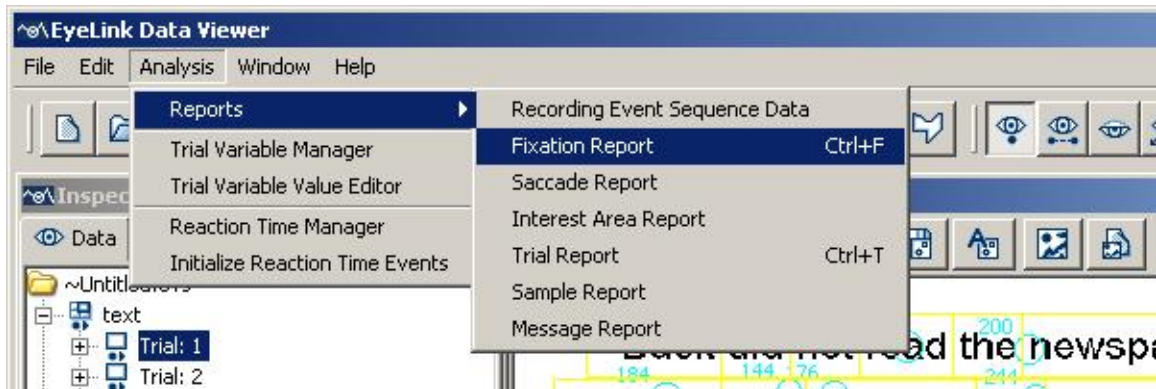


Figure 6-19. Creating a Fixation Report.

- 2) In the following output report dialog, Select the variables in the list of available output variables (left panel) and press ">>" to enter the variables into the list of selected output variables.

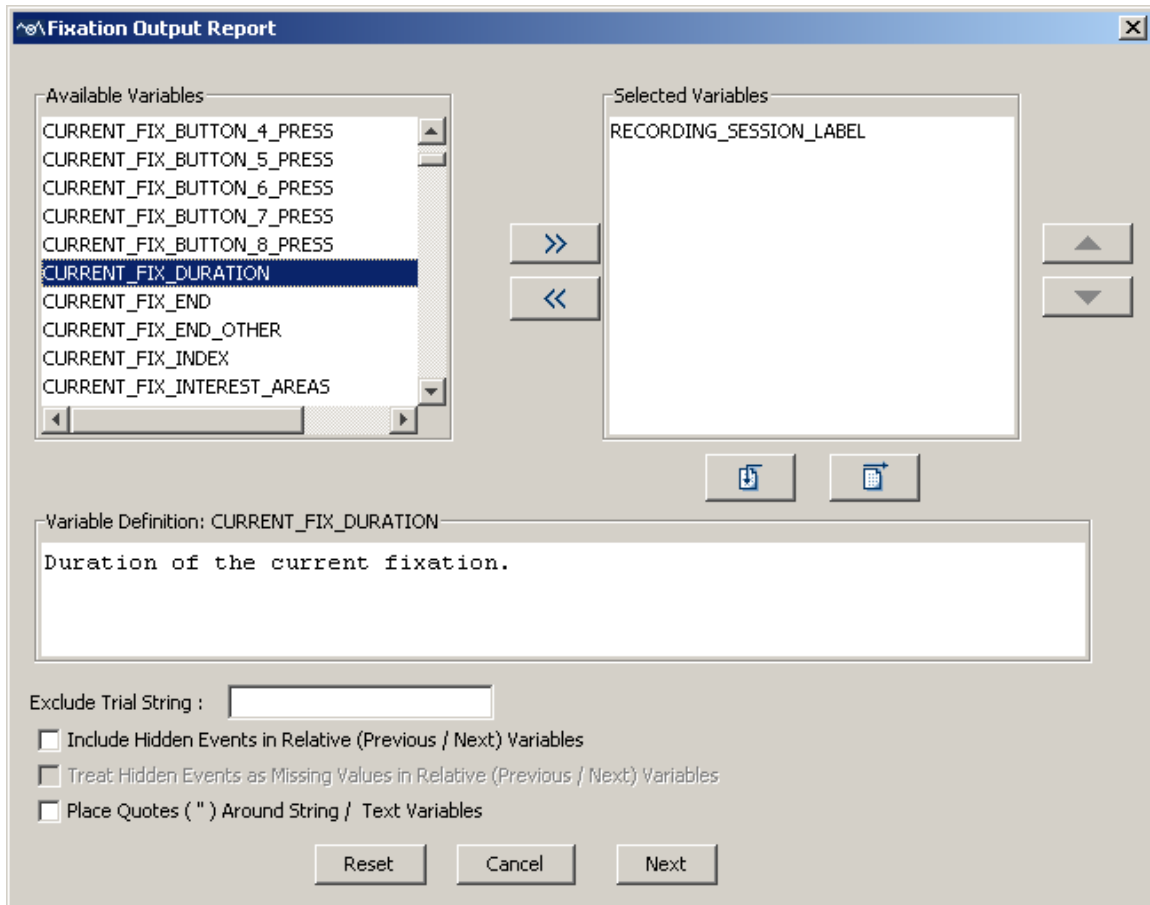




Figure 6-20. Variables in a Fixation Report.

- To remove variables from the output list, simply select the variables and press <<.
- To change the order in which the variables are listed, simply select one variable and use the ▲ or ▼ button to move it to the desired position. Repeat this step until all variables are in the right position.
- Data Viewer supports exporting and importing of variable selection across viewing sessions. For example, the user may go through each of the output reports to configure the selection of variables in an intended order. Export the variable selection into a property file by clicking on the "Export Variable Selection" button . For all future viewing sessions, the variable selection in the reports can be automated by importing the previously saved property file (by clicking on the "Import Variable Selection" button ).

3) Once all the desired variables have been entered into the output list, press NEXT button.

4) In the Save dialog box, browse to the directory in which you want to save the output report. The default output directory can be configured through "General Preferences".

5) In the File name text box, enter the name for the document and choose the appropriate file extension. Note, if the output report file extension set to "All Files" and the file name does not contain an extension, the file name is appended with a ".txt" by default.

6) Press the "Save" button.

6.6.2 Variables in the Fixation Report

Currently, the fixation report includes the following variables:

Field	Contents
RECORDING_SESSION_LABEL	Label of the data file
CURRENT_FIX_ADJUSTED	Whether the current fixation has been adjusted manually
CURRENT_FIX_BLINK_AROUND	Whether there is a blink preceding or following the current fixation. Possible values are "AFTER" (a blink follows the current fixation), "BEFORE" (a blink precedes the current fixation), "BOTH" (the current fixation is flanked by a preceding and a following blink), or "NONE" (no blink appears around the current fixation).
CURRENT_FIX_BUTTON_0_PRESS ... CURRENT_FIX_BUTTON_8_PRESS	Time from trial start if a specific button (0 - 8) has been pressed during or after fixation. If no button press is made from the start of the current fixation to the start of next fixation, missing value "." is assigned.
CURRENT_FIX_DURATION	Duration of the current fixation
* CURRENT_FIX_END	Trial time when the current fixation ends
* CURRENT_FIX_END_OTHER	Trial time when the current fixation ends (on the other eye in a binocular recording).
CURRENT_FIX_INDEX	Ordinal sequence of the current fixation in the trial
CURRENT_FIX_INTEREST_AREAS	Interest area in which the current fixation falls.
CURRENT_FIX_INTEREST_AREA_DWELL_TIME	Total amount of fixation dwell time (in msec) on the interest area to which the current fixation belongs.
CURRENT_FIX_INTEREST_AREA_FIX_COUNT	Total number of fixations on the interest area to which the current fixation belongs.
CURRENT_FIX_INTEREST_AREA_DATA	Reports the data (interest area type, dynamic/static interest area, and coordinate data) of the interest area to which the current fixation is assigned.

CURRENT_FIX_INTERESTAREA_GROUP	Group label for the interest area to which the current fixation is assigned.
CURRENT_FIX_INTERESTAREA_ID	The ID of the interest area in which the current fixation falls.
CURRENT_FIX_INTERESTAREA_INDEX	The index of the interest area in which the current fixation falls.
CURRENT_FIX_INTERESTAREA_LABEL	Label for the interest area to which the current fixation is assigned.
CURRENT_FIX_INTERESTAREA_PIXEL_AREA	Pixel area of the interest area to which the current fixation is assigned.
CURRENT_FIX_INTERESTAREA_RUN_ID	The ordinal sequence of the current run ² of fixation(s) made towards the current interest area. If the current interest area does not have a previous run of fixation(s) on it, this variable will have a value of 1.
CURRENT_FIX_INTERESTAREA_X_OFFSET	Horizontal offset of the current fixation relative to the center of the interest area. If the fixation falls into multiple interest areas, returns the offset value relative to the first relevant interest area only. Returns a missing value (.) for freehand interest areas.
CURRENT_FIX_INTERESTAREA_Y_OFFSET	Vertical offset of the current fixation relative to the center of the interest area.
CURRENT_FIX_IS_RT_END	Whether the current fixation is the end event of RT definition
CURRENT_FIX_LABEL	Label of the current fixation.
CURRENT_FIX_MSG_COUNT	Number of visible messages associated with the current fixation event ¹ . This number will not exceed the “Maximum Message Variables for Eye Event” set in the Output/Analysis Preference Settings.
CURRENT_FIX_MSG_TEXT_?	Text string for messages associated with the current fixation ¹ .
CURRENT_FIX_MSG_TIME_?	Trial time of the message ¹ .
CURRENT_FIX_NEAREST_INTEREST_AREA	Nearest interest area the current fixation is assigned to
CURRENT_FIX_NEAREST_INTEREST_AREA_DISTANCE	Distance (in degrees of visual angle) between the current fixation point and the center of the nearest interest area the current fixation is associated with
CURRENT_FIX_NEAREST_INTEREST_AREA_LABEL	Label for the nearest interest area the current fixation is associated with
CURRENT_FIX_PUPIL	Pupil size of the current fixation

CURRENT_FIX_REFIX_INTEREST_AREA	Whether this trial has a previous fixation in a different run ² that had the same interest area as the current fixation. If so, the CURRENT_FIX_INDEX value of the very first fixation in that interest area is given.
CURRENT_FIX_REFIX_PREV_INTEREST_AREA	Whether this trial has a previous fixation in an IA with higher IA ID than current fix IA ID (see run ² definition).
CURRENT_FIX_RUN_DWELL_TIME	Total amount of dwell time (in msec) in the current run ² of fixations.
CURRENT_FIX_RUN_INDEX	Ordinal sequence of the current fixation in the current run ² of fixations. The first fixation in this run has a RUN_INDEX of 1 and the second one has RUN_INDEX value of 2, and so on.
CURRENT_FIX_RUN_SIZE	Number of fixations within the current run ² .
* CURRENT_FIX_START	Trial time when the current fixation starts
* CURRENT_FIX_START_OTHER	Trial time when the current fixation starts (on the other eye in a binocular recording).
CURRENT_FIX_TRIAL_SPAN	Whether the fixation starts before the trial starts and ends after the trial starts, or starts before the trial ends and ends after the trial ends
CURRENT_FIX_X	X coordinate of the current fixation
CURRENT_FIX_X_OTHER	X coordinate of the current fixation (on the other eye in a binocular recording)
CURRENT_FIX_X_RESOLUTION	Horizontal angular resolution (in screen pixels per degree) for the current fixation.
CURRENT_FIX_Y	Y coordinate of the current fixation
CURRENT_FIX_Y_OTHER	Y coordinate of the current fixation (on the other eye in a binocular recording).
CURRENT_FIX_Y_RESOLUTION	Vertical angular resolution (in screen pixels per degree) for the current fixation.
	Note: Similar analyses are done for the previous fixation (variables beginning with PREVIOUS_FIX) and for the next fixation (variables beginning with NEXT_FIX).
DATA_FILE	File name of the recording;
EYE_USED	Which eye's data (LEFT or RIGHT) is used to create this report?
IP_END_TIME	End time (in milliseconds since EyeLink tracker was activated) of the interest period.
IP_LABEL	Label of the current interest period selected in the viewing session.

IP_START_TIME	Start time (in milliseconds since EyeLink tracker was activated) of the interest period.
LAST_BUTTON_PRESSED	ID of the pressed button
* LAST_BUTTON_PRESS_TIME	Trial time (relative to "TRIAL_START_TIME") when the last button is pressed
LAST_BUTTON_RELEASED	ID of the released button
*LAST_BUTTON_PRELEASE_TIME	Trial time (relative to "TRIAL_START_TIME") when the last button is released
* LAST_BUTTON_TIME	Trial time (relative to "TRIAL_START_TIME") when the last button is pressed or released.
NEXT_FIX_ANGLE, PREVIOUS_FIX_ANGLE	Angle between the horizontal plane and the line connecting the current fixation and the previous/next fixation (see 5-2)
NEXT_FIX_DIRECTION, PREVIOUS_FIX_DIRECTION	Direction (Left, Right, Top, Bottom), relative to the current fixation, in which the next/previous fixation is located
NEXT_FIX_DISTANCE, PREVIOUS_FIX_DISTANCE	Distance (in degrees of visual angle) between the current fixation and the next/previous fixation in visual degrees
NEXT_SAC_AMPLITUDE	Amplitude of the following saccade in degrees of visual angle
NEXT_SAC_ANGLE	Angle between the horizontal plane and the direction of the next saccade (see 5-2)
NEXT_SAC_AVG_VELOCITY	Average velocity of the next saccade
NEXT_SAC_CONTAINS_BLINK	Whether the next saccade contains blink(s).
NEXT_SAC_DIRECTION	Direction (Left, Right, Top, Bottom), relative to the current fixation, in which the next saccade is aiming
NEXT_SAC_DURATION	Duration of the next saccade in milliseconds
NEXT_SAC_END_INTEREST AREAS	Interest area in which the end point of the next saccade falls
NEXT_SAC_END_INTEREST_A REA_ID	The ID of the interest area to which the end point of the next saccade is assigned.
NEXT_SAC_END_INTEREST_A REA_INDEX	Interest area to which the end point of the next saccade is assigned.
NEXT_SAC_END_INTEREST_A REA_LABEL	Label for the interest area to which the end position of the next saccade is assigned.
* NEXT_SAC_END_TIME	Trial time when the next saccade ends
NEXT_SAC_END_X	X coordinates of the end point for the next saccade
NEXT_SAC_END_X_RESOLUTION	Horizontal angular resolution (in screen pixels per degree) at the end of the saccade. If samples are loaded into the viewing session, this takes the horizontal resolution of the last sample in the

	saccade; otherwise, the horizontal resolution of the next fixation is used.
NEXT_SAC_END_Y	Y coordinates of the end point for the next saccade
NEXT_SAC_END_Y_RESOLUTION	Vertical angular resolution (in screen pixels per degree) at the end of the saccade. If samples are loaded into the viewing session, this takes the vertical resolution of the last sample in the saccade; otherwise, the vertical resolution of the next fixation is used.
NEXT_SAC_INDEX	Ordinal index of the next saccade
NEXT_SAC_IS_RT_END	Whether the next saccade is the end event of the RT definition
NEXT_SAC_LABEL	Label of the next saccade.
NEXT_SAC_MSG_COUNT	Number of visible messages associated with the next saccade event ¹ . This number will not exceed the "Maximum Message Variables for Eye Event" set in the Output/Analysis Preference Settings.
NEXT_SAC_MSG_TEXT	Text string of the messages associated with the next saccade ¹ .
NEXT_SAC_MSG_TIME	Trial time ¹ of the message ¹ .
NEXT_SAC_NEAREST_END_INTEREST_AREA	Nearest interest area to which the end point of the next saccade is assigned
NEXT_SAC_NEAREST_END_INTEREST_AREA_LABEL	Label for the nearest interest area to which the end point of the next saccade is assigned
NEXT_SAC_NEAREST_START_INTEREST_AREA	Nearest interest area to which the start point of the next saccade is assigned
NEXT_SAC_NEAREST_START_INTEREST_AREA_LABEL	Label for the nearest interest area to which the start point of the next saccade is assigned
NEXT_SAC_PEAK_VELOCITY	Peak values of gaze velocity (in visual degrees per second) of the next saccade
NEXT_SAC_START_INTEREST_AREAS	Interest area in which the start point of the next saccade falls
NEXT_SAC_START_INTEREST_AREA_ID	The ID of the interest area to which the start point of the next saccade is assigned.
NEXT_SAC_START_INTEREST_AREA_INDEX	Interest area to which the start point of the next saccade is assigned.
NEXT_SAC_START_INTEREST_AREA_LABEL	Label for the interest area to which the start position of the next saccade is assigned.
* NEXT_SAC_START_TIME	Trial time when the next saccade ends
NEXT_SAC_START_X	X coordinates of the start point for the next saccade
NEXT_SAC_START_X_RESOLUTION	Horizontal angular resolution (in screen pixels per

TION	degree) at the start of the saccade. If samples are loaded into the viewing session, this takes the horizontal resolution of the first sample in the saccade; otherwise, the horizontal resolution of the current fixation is used.
NEXT_SAC_START_Y	Y coordinates of the start point for the next saccade
NEXT_SAC_START_Y_RESOLUTION	Vertical angular resolution (in screen pixels per degree) at the start of the saccade. If samples are loaded into the viewing session, this takes the vertical resolution of the first sample in the saccade; otherwise, the vertical resolution of the current fixation is used.
	Note: Similar analyses are done for the previous saccade (variables beginning with PREVIOUS_SAC)
TRIAL_FIXATION_TOTAL	Total number of fixations in a trial
TRIAL_INDEX	Sequential order of the trial in the recording.
TRIAL_LABEL	Label of the trial
* TRIAL_START_TIME	The start time (in milliseconds since EyeLink tracker was activated) of the trial.
VIDEO_FRAME_INDEX_START	Index of the video frame that was visible at the start of the eye event. Note that all “VIDEO_” variables are applicable to SceneLink recordings or some Experiment Builder recordings only. For SceneLink recordings, please check out the “Data Loading” Preference Settings.
VIDEO_FRAME_INDEX_END	Index of the video frame that was visible at the end of the eye event.
VIDEO_NAME_START	Name of the video frame that was visible at the start of the eye event.
VIDEO_NAME_END	Name of the video frame that was visible at the end of the eye event.

In addition to these default variables, other user-defined variables are also listed if they have been defined in the EDF file (see section 7.2.1 on creating trial variable labels) or created during the viewing session (see section 6.1.1 on trial variable manager).

Please note that the data output in the fixation report depends on the setting of interest period filter (see section 6.4.2 or Figure 6-18) – only those fixation and saccade events falling within the active interest period will be reported. If the "Trimmed Spanned Fixation Duration" setting of the "Data Filter" preference is enabled, the duration, start

time, and end time of the fixations that overlap with the start or end of the interest period will be adjusted accordingly. If a valid reaction time definition is found in a trial, the TRIAL_START_TIME will be set as the start as the reaction time definition; otherwise this will be set as the start of trial recording (see section 6.5.3 on discussion of three time periods used in Data Viewer). As a result, the values of CURRENT_FIX_END and CURRENT_FIX_START, which are calculated relative to TRIAL_START_TIME, will be different depending on whether a valid reaction time definition is found for the trial.

Note ¹: Messages written in the EDF file will be associated with an event closest in time. So, for a fixation report, the message could appear during a preceding saccade (PREVIOUS_SAC_MSG_*) or a following saccade (NEXT_SAC_MSG_*). In very rare cases, the message can be written to both saccade and fixation events if the message was written out after the last sample of a fixation (or saccade) but before the first sample of the following saccade (or fixation).

Note ²: A run of fixations refers to a group of consecutive fixations that are directed towards the same interest area. Thus, for example, the interest areas of 10 fixations are: A A B A A A C A E D. The run count will be 7. The following table summarizes the values for some relevant variables.

Variables	A	A	B	A	A	A	C	A	E	D
Run (not listed in the variable list)	1	1	2	3	3	3	4	5	6	7
CURRENT_FIX_INDEX	1	2	3	4	5	6	7	8	9	10
CURRENT_FIX_INTEREST_A REA_INDEX	1	1	2	1	1	1	3	1	5	4
CURRENT_FIX_INTEREST_A REA_FIX_COUNT	6	6	1	6	6	6	1	6	1	1
CURRENT_FIX_INTEREST_A REA_RUN_ID	1	1	1	2	2	2	1	3	1	1
CURRENT_FIX_REFIX_INTER EST_AREA	0	0	0	1	1	1	0	1	0	0
CURRENT_FIX_REFIX_PREV_ INTEREST_AREA	0	0	0	1	1	1	0	1	0	1
CURRENT_FIX_RUN_INDEX	1	2	1	1	2	3	1	1	1	1
CURRENT_FIX_RUN_SIZE	2	2	1	3	3	3	1	1	1	1

6.6.3 Configuring the Fixation Report

In addition to excluding some trials by entering a string in the “Exclude Trial String” edit box, the fixation report (similarly the saccade report and the interest area report) also

allows the inclusion of hidden events in the relative event variables [PREVIOUS_ and NEXT_]. This can be done by clicking the “Include Hidden Events in Relative Variables” check box. If the hidden events are included in the analysis, the user can further decide whether or not to treat these events as missing data.

6.7 Saccade Report

Saccade report creates a columnar output of saccade events in the data file. Each row of the report represents a saccade event. Depending on the variables selected for output, each row can contain information on the previous fixation and/or saccade, as well as the next fixation and/or saccade.

6.7.1 Obtaining a Saccade Report

To obtain a saccade report, adopt the same procedure as in obtaining the trial report, except from the menus, choose:

Analysis → Report → Saccade Report

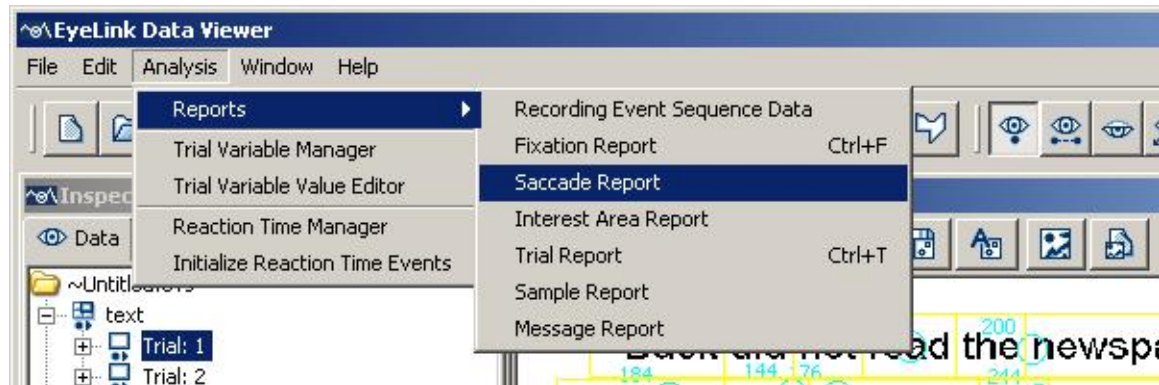


Figure 6-21. Creating a Saccade Report.

6.7.2 Variables in the Saccade Report

Currently, the saccade report includes the following saccade-related information: index, start time, end time, start *xy* coordinates, end *xy* coordinates, duration, amplitude, angle, direction, average velocity, peak velocity, interest areas of the current, previous, and next saccades, and whether these saccades have been selected for analysis. It also contains the following information on the previous fixation and the next fixation: angle, blink around, direction, distance, duration, start time, end time, pupil size, trial span, the interest area in which the fixation is located, the nearest interest area, and whether the fixation has been selected for analysis. In addition, the report contains labels of the data file, data file, and individual trials, as well as other user-defined variables (see section 6.6.2 Variables in the Fixation Report for the complete list of variables and their meanings).

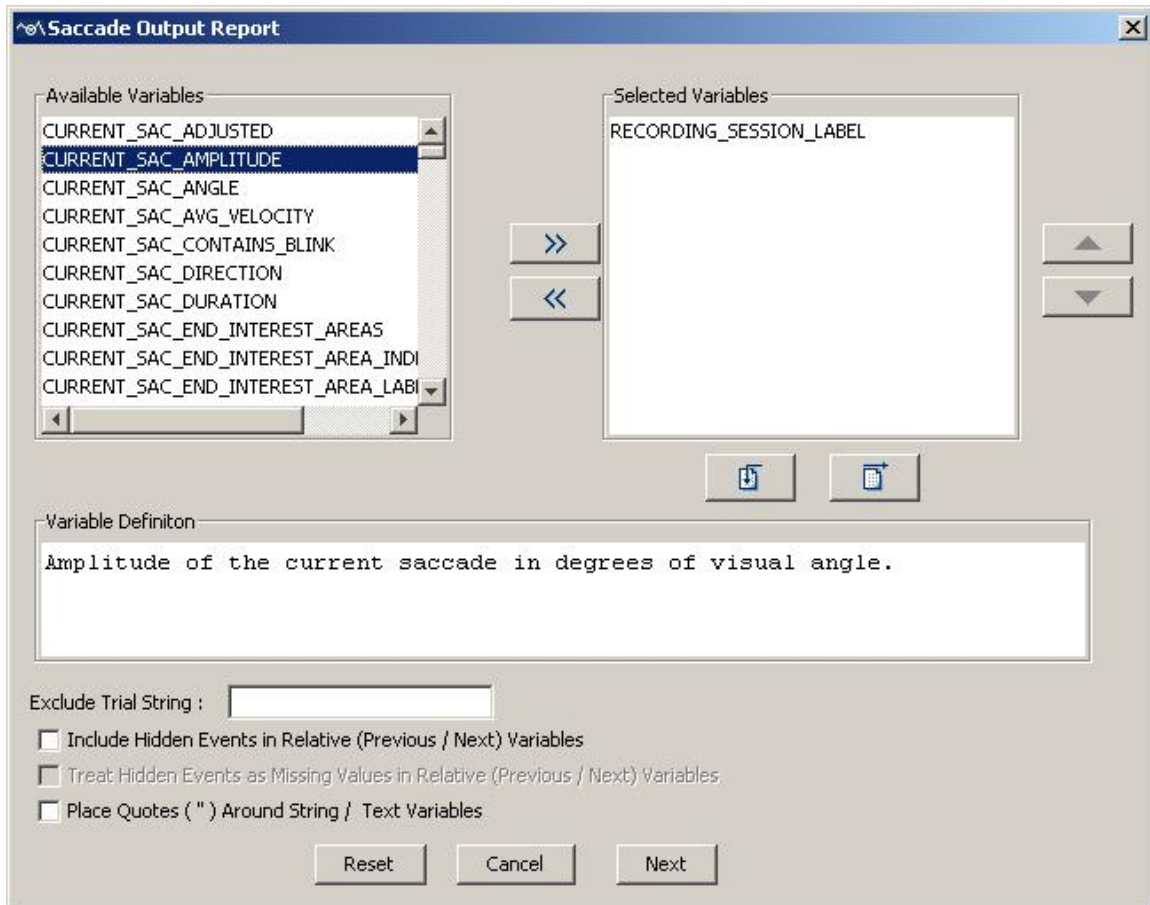


Figure 6-22. Variables in a Saccade Report.

6.8 Interest Area Report

Interest area report provides a columnar output of eye movement data for each interest area in a trial. Each row of the report represents an interest area, arranged in an ascending order of the interest area ID.

6.8.1 Obtaining an Interest Area Report

To obtain an interest area report,

- 1) From the application menus, choose:
Analysis -> Report -> Interest Area Report

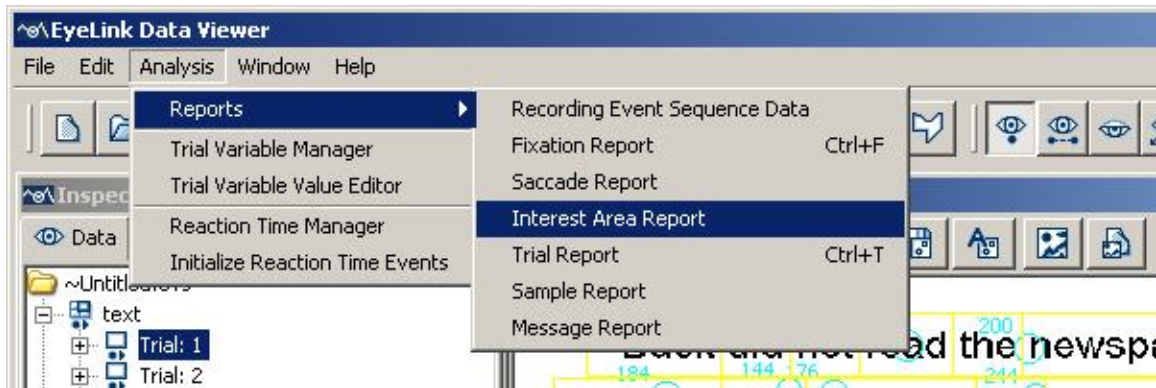


Figure 6-23. Creating an Interest Area Report.

2) In the following output report dialog, Select the variables in the list of available output variables (left panel) and press ">>" to enter the variables into the list of selected output variables.

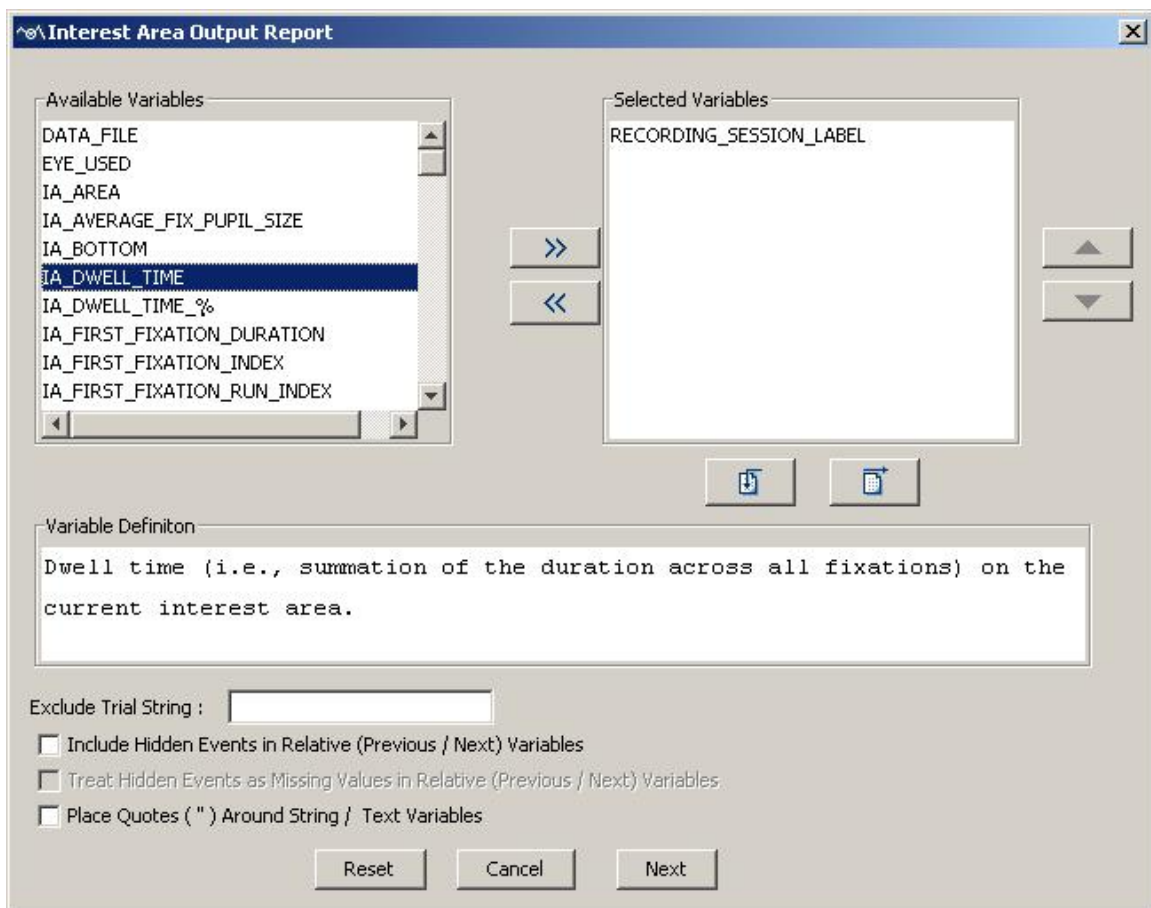




Figure 6-24. Variables in an Interest Area Report.

- To remove variables from the output list, simply select the variables and press <<.
- To change the order in which the variables are listed, simply select one variable and use the ▲ or ▼ button to move it to the desired position. Repeat this step until all variables are in the right position.
- Data Viewer supports exporting and importing of variable selection across viewing sessions. For example, the user may go through each of the output reports to configure the selection of variables in an intended order. Export the variable selection into a property file by clicking on the "Export Variable Selection" button . For all future viewing sessions, the variable selection in the reports can be automated by importing the previously saved property file (by clicking on the "Import Variable Selection" button ).

3) Once all the desired variables have been entered into the output list, press NEXT button.

4) In the Save dialog box, browse to the directory in which you want to save the output report. The default output directory can be configured through "General Preferences".

5) In the File name text box, enter the name for the document and choose the appropriate file extension. Note, if the output report file extension set to "All Files" and the file name does not contain an extension, the file name is appended with a ".txt" by default.

6) Press the "Save" button.

Data Viewer allows the user to configure the interest area report. The user can exclude some trials by entering a string in the "Exclude Trial String" edit box. This string should be one message or part of a message within the scope of the trial recording that uniquely defines that subset of trials to be excluded. In addition, to enable correct loading of string/text variables in some statistics software, the Viewer also has an option of adding a pair of quotation marks for such variables.

6.8.2 Variables in the Interest Area Report

Currently, the interest area report provides information for the following variables:

Field	Contents
RECORDING_SESSION_LABEL	Label of the data file
DATA_FILE	File name of the recording
EYE_USED	Which eye's data (LEFT or RIGHT) is used to create this report?
IA_AREA	Pixel area for the current interest area.
IA_AVERAGE_FIX_PUPIL_SIZE	Average pupil size across all fixations in the interest area.

IA_BOTTOM	Bottom side pixel position of non-freehand IA.
IA_DWELL_TIME	Dwell time (i.e., summation of the duration across all fixations) on the current interest area
IA_DWELL_TIME_%	Percentage of trial time spent on the current interest area
IA_DYNAMIC	Whether the current interest area is a dynamic one (true or false).
IA_END_TIME	Time (relative to the start of the current interest period) when the last instance of the current interest area ends. This variable is applicable to dynamic interest area only.
IA_FIRST_FIXATION_DURATION	Duration of the first fixation event that was within the current interest area.
IA_FIRST_FIXATION_INDEX	Ordinal sequence of the first fixation that was within the current interest area.
IA_FIRST_FIXATION_RUN_INDEX	This counts how many runs of fixations have occurred when a first fixation is made to an interest area. The current run is also included in the tally.
* IA_FIRST_FIXATION_TIME	Start time of the first fixation to enter the current interest area.
IA_FIRST_FIXATION_X	The X position of the first fixation event that was within the current interest area.
IA_FIRST_FIXATION_Y	The Y position of the first fixation event that was within the current interest area.
IA_FIRST_FIX_PROGRESSIVE	Checks whether later interest areas have been visited before the first fixation enters the current interest area. 1 if NO higher IA ID in earlier fixations before the first fixation in the current interest area; 0 otherwise. This measure is useful in reading to check whether the first run of fixations in this interest area is in fact first-pass fixations.
IA_FIRST_RUN_DWELL_TIME	Dwell time (i.e., summation of the duration across all fixations) of the first run within the current interest area.
* IA_FIRST_RUN_END_TIME	End time of the first run of fixations in the current interest area.
IA_FIRST_RUN_FIXATION_%	Percentage of all fixations in a trial falling in the first run of the current interest area.
IA_FIRST_RUN_FIXATION_COUNT	Number of all fixations in a trial falling in the first run of the current interest area.
* IA_FIRST_RUN_START_TIME	Start time of the first run of fixations in the current interest area.

IA_FIRST_SACCADE_AMPLITUDE	Amplitude (in degree of visual angle) of the first saccade entering into the current interest area.
IA_FIRST_SACCADE_ANGLE	Angle between the horizontal plane and the direction of the first saccade entering into the current interest area.
IA_FIRST_SACCADE_INDEX	Ordinal index of the saccade in a trial which first landed within the current interest area.
*IA_FIRST_SACCADE_END_TIME	End time of the saccade that first landed within the current interest area.
*IA_FIRST_SACCADE_START_TIME	Start time of the saccade that first landed within the current interest area.
IA_FIXATION_%	Percentage of all fixations in a trial falling in the current interest area
IA_FIXATION_COUNT	Total fixations falling in the interest area
IA_FSA_COUNT	The number of fixations (fixation N) which started in the current row of interest area, with fixation N + fixation_skip_count ending in the current column of interest area.
IA_FSA_DURATION	The summed duration for all fixations (fixation N) which started in the current row of interest area, with fixation N + fixation_skip_count ending in the current column of interest area.
IA_GROUP	Group label for the current interest area.
IA_ID	Ordinal ID of the current interest area
IA_INSTANCE_COUNT	The total number of individual instances of the current dynamic interest area.
IA_LABEL	Label for the current interest area
IA_LAST_FIXATION_DURATION	Duration of the last fixation event that was within the current interest area.
IA_LAST_FIXATION_RUN	Index of run that the last fixation is in.
*IA_LAST_FIXATION_TIME	Start time of the last fixation to enter the current interest area.
IA_LAST_FIXATION_X	The X position of the last fixation event that was within the current interest area.
IA_LAST_FIXATION_Y	The Y position of the last fixation event that was within the current interest area.
IA_LAST_RUN_DWELL_TIME	Dwell time (i.e., summation of the duration across all fixations) of the last run of fixations within the current interest area.
IA_LAST_RUN_END_TIME	End time of the last run of fixations in the current interest area.
IA_LAST_RUN_FIXATION_%	Percentage of all fixations in a trial falling in the

	last run in the current interest area.
IA_LAST_RUN_FIXATION_COUNT	Number of all fixations in a trial falling in the last run in the current interest area.
IA_LAST_RUN_START_TIME	Start time of the last run of fixations in the current interest area.
IA_LAST_SACCADE_AMPLITUDE	Amplitude (in degree of visual angle) of the last saccade entering into the current interest area.
IA_LAST_SACCADE_ANGLE	Angle between the horizontal plane and the direction of the last saccade entering into the current interest area.
*IA_LAST_SACCADE_END_TIME	End time of the saccade that last landed within the current interest area.
IA_LAST_SACCADE_INDEX	Ordinal index of the saccade in a trial which last landed within the current interest area.
*IA_LAST_SACCADE_START_TIME	Start time of the saccade that last landed within the current interest area.
IA_LEFT	Left side pixel position of non-freehand IA.
IA_LEGAL	A trial is considered legal (i.e., IA_LEGAL = 1) only if the first fixation in 'interest area N' was preceded by a fixation in interest area (N-1). This variable is relevant for experiments that run in a contingent display paradigm, as it ensures that readers could see the target area before actually landing in it.
IA_MAX_FIX_PUPIL_SIZE	Maximum pupil size among all fixations in the interest area. Note this measure does not represent the maximum pupil size within a fixation.
IA_MIN_FIX_PUPIL_SIZE	Minimum pupil size among all fixations in the interest area. Note this measure does not represent the minimum pupil size within a fixation.
IA_POINTS	List of x,y points for freehand Interest Areas.
IA_REGRESSION_IN	Whether the current interest area received at least one regression from later interest areas (e.g., later parts of the sentence). 1 if interest area was entered from a higher IA_ID (from the right in English); 0 if not.
IA_REGRESSION_IN_COUNT	Number of times interest area was entered from a higher IA_ID (from the right in English).
IA_REGRESSION_OUT	Whether regression(s) was made from the current interest area to earlier interest areas (e.g., previous parts of the sentence) prior to leaving that interest area in a forward direction. 1 if a saccade exits the

	current interest area to a lower IA_ID (to the left in English) before a later interest area was fixated; 0 if not.
IA_REGRESSION_OUT_FULL	Whether regression(s) was made from the current interest area to earlier interest areas (e.g., previous parts of the sentence). 1 if a saccade exits the current interest area to a lower IA_ID (to the left in English); 0 if not. Note that IA_REGRESSION_OUT only considers first-pass regressions whereas IA_REGRESSION_OUT_FULL considers all regressions, regardless whether later interest areas have been visited or not.
IA_REGRESSION_OUT_FULL_COUNT	Number of times interest area was exited to a lower IA_ID (to the left in English).
IA_REGRESSION_OUT_COUNT	Number of times interest area was exited to a lower IA_ID (to the left in English) before a higher IA_ID was fixated in the trial.
IA_REGRESSION_PATH_DURATION	The summed fixation duration from when the current interest area is first fixated until the eyes enter an interest area with a higher IA_ID.
IA_RIGHT	Right side pixel position of non-freehand IA.
IA_REGRESSION_PATH_DURATION	The summed fixation duration from when the current interest area is first fixated until the eyes enter an interest area with a higher IA_ID.
IA_RUN_COUNT	Number of times the Interest Area was entered and left (runs)
IA_SECOND_FIXATION_DURATION	Duration of the second fixation in IA, regardless of run.
IA_SECOND_FIXATION_RUN	Index of run that the second fixation is in.
* IA_SECOND_FIXATION_TIME	Time of the second fixation in IA, regardless of run.
IA_SECOND_FIXATION_X	The X position of the second fixation event that was within the current interest area.
IA_SECOND_FIXATION_Y	The Y position of the second fixation event that was within the current interest area.
IA_SECOND_RUN_DWELL_TIME	Dwell time (i.e., summation of the duration across all fixations) of the second run of fixations within the current interest area.
* IA_SECOND_RUN_END_TIME	End time of the second run of fixations in the current interest area.
IA_SECOND_RUN_FIXATION	Percentage of all fixations in a trial falling in the

%	second run of the current interest area.
IA_SECOND_RUN_FIXATION_COUNT	Number of all fixations in a trial falling in the second run in the current interest area.
*IA_SECOND_RUN_START_TIME	Start time of the second run of fixations in the current interest area.
IA_SELECTIVE_REGRESSION_PATH_DURATION	Duration of fixations and refixations of the current interest area before the eyes enter an interest area with a higher ID.
IA_SKIP	An interest area is considered skipped (i.e., IA_SKIP = 1) if no fixation occurred in first-pass reading.
IA_SPILLOVER	The duration of the first fixation made on 'interest area (N+1)' after leaving 'interest area N' in first pass.
IA_START_TIME	Time (relative to the start of the current interest period) when the first instance of the current interest area starts. This variable is applicable to dynamic interest area only.
IA_THIRD_FIXATION_DURATION	Duration of the third fixation in IA, regardless of run.
IA_THIRD_FIXATION_RUN	Index of run that the third fixation is in.
*IA_THIRD_FIXATION_TIME	Time of the third fixation in IA, regardless of run.
IA_THIRD_FIXATION_X	The X position of the third fixation event that was within the current interest area.
IA_THIRD_FIXATION_Y	The Y position of the third fixation event that was within the current interest area.
IA_THIRD_RUN_DWELL_TIME	Dwell time (i.e., summation of the duration across all fixations) of the third run of fixations within the current interest area.
*IA_THIRD_RUN_END_TIME	End time of the third run of fixations in the current interest area.
IA_THIRD_RUN_FIXATION_COUNT	Number of all fixations in a trial falling in the third run in the current interest area.
IA_THIRD_RUN_FIXATION_%	Percentage of all fixations in a trial falling in the third run of the current interest area.
*IA_THIRD_RUN_START_TIME	Start time of the third run of fixations in the current interest area.
IA_TYPE	Type of interest area (rectangular, elliptical, or freehand)
IA_TOP	Top side pixel position of non-freehand IA
IP_END_TIME	End time (in milliseconds since EyeLink tracker

	was activated) of the interest period.
IP_LABEL	Label of the current interest period selected in the viewing session.
IP_START_TIME	Start time (in milliseconds since EyeLink tracker was activated) of the interest period.
REPORTING_METHOD	Whether a fixation-based or sample-based calculation is used in creating the interest area report. The reporting method can be configured through "Output/Analysis Preference -> Perform Sample-based Calculation".
TRIAL_DWELL_TIME	Dwell time (i.e., summation of all fixation durations) for the whole trial
TRIAL_FIXATION_COUNT	Total number of fixations in the trial
TRIAL_IA_COUNT	Reports the total number of interest areas found in the trial.
TRIAL_INDEX	Sequential order of the trial in the recording.
TRIAL_LABEL	Label of the trial
* TRIAL_START_TIME	The start time (in milliseconds since EyeLink tracker was activated) of the trial.
TRIAL_TOTAL_VISITED_IA_COUNT	Reports total number of unique interest areas visited in the trial.

In addition, other user-defined trial variable labels will also be listed. Note: since version 1.3 of Data Viewer, all eye event start and end times in the interest area report are made relative to trial start instead. The absolute time of eye events can be obtained by adding "TRIAL_START_TIME".

6.8.3 Using Interest Area Report

The data output in the interest area report depends on the setting of the interest period filter (see section 6.4.2 or Figure 6-18) – only those fixation and saccade events falling within the active interest period will be reported. If a valid reaction time definition is found in a trial, the TRIAL_START_TIME will be set as the start as the reaction time definition; otherwise this will be set as the start of trial recording (see section 6.5.3 on discussion of three time periods used in Data Viewer). As a result, the values of all of variables marked with a "*" symbol, which are calculated relative to TRIAL_START_TIME, will be different depending on whether a valid reaction time is found in the trial. If the "Trimmed Spanned Fixation Duration" setting of the "Data Filter" preference is enabled, the duration, start time, and end time of the fixations that overlap with the start or end of the interest period will be adjusted accordingly.

Some variables in the interest area report may be influenced by the outlier fixations that do not belong to any interest area. For example, the presence of an outlier fixation may break the run of fixations on a particular interest area and therefore shortens the measure of IA_FIRST_RUN_DWELL_TIME and increases the IA_RUN_COUNT value. Therefore, users should ensure that all fixations are assigned to a particular interest area before creating an interest area report, if "run"-related measures are important. Users can either manually move the position of those outlier fixations or, more conveniently, enable the "Use Nearest Interest Area for Outlier Fixations" option of the Output/Analysis preference settings.

Interest area report was designed to support analysis of data from different research paradigms and, therefore, the user only needs to use a small portion of variables in one output. **If you have lots of interest areas in a trial, including IA_FSA_COUNT and IA_FSA_DURATION in the output report may take a very long time for the Viewer to generate the report.** A progress dialog (see Figure 6-25) will be displayed while creating the report. The user may abort the process by pressing the "Cancel" button.


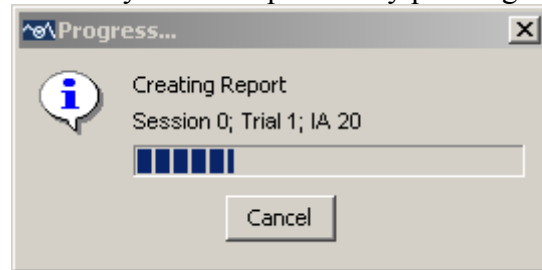


Figure 6-25. Progress Dialog Box Displayed when Creating the Interest Area Report

6.8.3.1 Fixation Sequence Analysis

Variables IA_FSA_COUNT and IA_FSA_DURATION are used primarily for fixation sequence analysis (i. e., examination of the frequency of saccades directed from the current interest area to all interest areas). If there are N interest areas in a trial, this will generate N columns in the output report, with column labels being IA_FSA_COUNT_1 to IA_FSA_COUNT_N for counts and IA_FSA_DURATION_1 to IA_FSA_DURATION_N for duration outputs. Each column in the output file reports number of fixations coming to one particular interest area from each of all possible interest areas. Each row reports number of fixations starting from the current interest area and ending at each of the interest areas (when "Fixation Skip Count" in the Output/Analysis preference settings is set to 1).

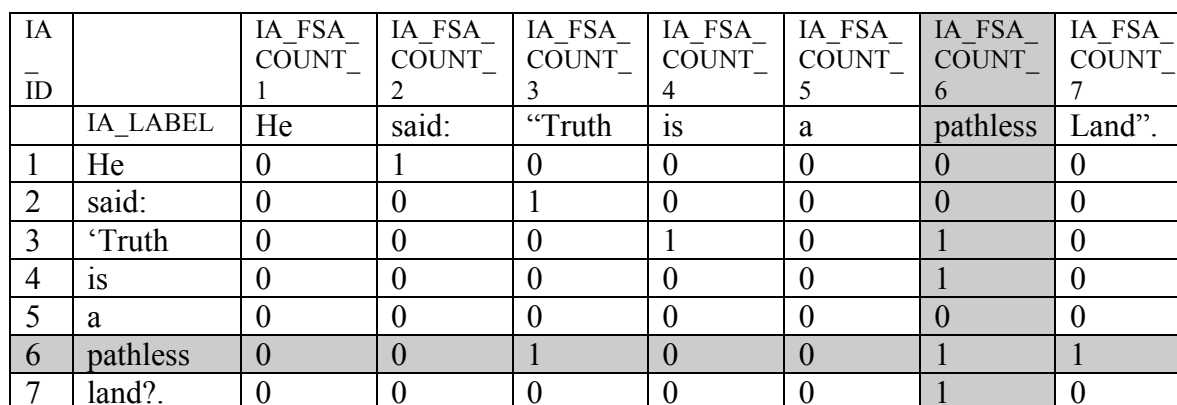

Please note that the user should also check the setting for “Fixation Skip Count” in the Output/Analysis preference settings. The fixation skip count defines the number (should be an integer no less than 1) of fixations to skip when looking for the next interest area to use in the fixation sequence analysis variables of the Interest Area report. The default setting is 1 and uses the next available fixation for the sequence analysis (from the current interest area to all interest area). A value of 2 uses the fixation following the next one for the analysis, and so on.

A lot of the variables in the interest area report are intended for the regression analysis, especially in the reading research. The following example illustrates the use the interest area report in the regression analysis.

IA_ID	1	2	3	4	5	6	7
IA_LABEL	He	said:	“Truth	is	a	pathless	Land”.
IA_DWELL_TIME	F1	F2	F3 + F6	F7	0	F4 + F5 + F8 + F10	F9
IA_FIXATION COUNT	1	1	2	1	0	4	1
IA_RUN_COUNT	1	1	2	1	.	3	1
IA_FIRST_FIXATION INDEX	1	2	3	7	.	4	9
IA_FIRST_FIXATION _RUN_INDEX	1 [1]	2 [1, 2]	3 [1, 2, 3]	6 [1, 2, 3, 6, 3, 4]	.	4 [1, 2, 3, 6]	8 [1, 2, 3, 6, 3, 4, 6, 7]
IA_FIRST_FIXATION _VISITED_IA_COUNT	0 []	1 [1]	2 [1, 2]	4 [1, 2, 3, 6]	.	3 [1, 2, 3]	5 [1, 2, 3, 6, 4]
IA_FIRST_FIX PROGRESSIVE	1	1	1	0	.	1	1
IA_FIRST_FIXATION DURATION	F1	F2	F3	F7	.	F4	F9
IA_FIRST_RUN DWELL_TIME	F1	F2	F3	F7	.	F4 + F5	F9
IA_SECOND_FIXATION DURATION	.	.	F6	.	.	F5	.
IA_SECOND_FIXATION RUN	.	.	2	.	.	1	.
IA_SECOND_RUN DWELL_TIME	.	.	F6	.	.	F8	.
IA_REGRESSION_IN	0	0	1	0	.	1	0
IA_REGRESSION_IN COUNT	0	0	1	0	.	1	0
IA_REGRESSION OUT	0	0	0	0	.	1	1
IA_REGRESSION OUT_COUNT	0	0	0	0	.	1	1
IA_REGRESSION _PATH_DURATION	F1	F2	F3	F7	.	F4 + F5 + F6 + F7 + F8	F9
IA_SELECTIVE _REGRESSION_PATH DURATION	F1	F2	F3	F7	.	F4 + F5 + F8	F9

Take interest area #6 (“pathless”) for example. IA_REGRESSION_PATH_DURATION includes the first pass fixations on the interest area (F4 + F5), time spent in previous parts of the sentence following regressive eye movements (F6 + F7), and time due to

refixations coming from the left before the eyes move past the interest area (F8). The calculation of `IA_SELECTIVE_REGRESSION_PATH_DURATION` only includes the first run fixation on the interest area (F4+F5) and the refixations on the interest area (F8) before the eyes enter next interest area with a higher `IA_ID` ("land").

`IA_FIRST_FIX_PROGRESSIVE` checks whether later interest areas have been visited before the first fixation enters the current interest area. Therefore, this variable would be 0 for the interest area "is" because it was fixated on Fixation 7 while Fixations 4 and 5 were on later part of the sentence ("pathless"). `IA_REGRESSION_OUT` checks whether at least one regression was made from the current interest area to previous parts of the sentence prior to leaving that interest area in a forward direction. Therefore, the transition from F5 → F6 makes this flag true for the interest area #6 (labeled as "pathless"). `IA_REGRESSION_IN` checks whether a given interest area received at least one regression from later parts of the sentence. Therefore, this is true for both interest areas #3 "truth" (F5 → F6) and #6 "pathless" (F9 → F10).

6.9 Sample Output Report

Sample output report provides a columnar output of eye movement data for each eye sample in a trial. The sample report option will not be available if the samples are not loaded into a viewing session. As in the other reports, sample output report only outputs samples that fall within the active interest period for the trial.

6.9.1 Loading Samples into a Viewing Session

Version 1.7 of Data Viewer introduced a change to optimize the viewer performance by not loading the samples by default (see revision history). To make sample data available to a viewing session, you will need to first change the data loading preferences and then reload the EDF file.

- 1) Before loading any EDF files, please go to "Start -> Programs -> SR Research -> EyeLink -> Data Viewer" to start the software.
- 2) Once the application starts up, click "Preferences" tab and go to "Data Loading" preferences. Check the "Load Samples" option. If you want this changes to be persistent over future viewing sessions, you may consider saving this as default setting. This can be done by selecting the topmost preference tree node, clicking the right mouse buttons and selecting "Save Properties as Defaults".
- 3) Once you have done step 2), this will make the properties persist over future data viewing sessions. Now, start loading in the EDF files.

6.9.2 Obtaining a Sample Output Report

To obtain an output report for samples,

- 1) From the application menus, choose:
Analysis -> Report -> Sample Report

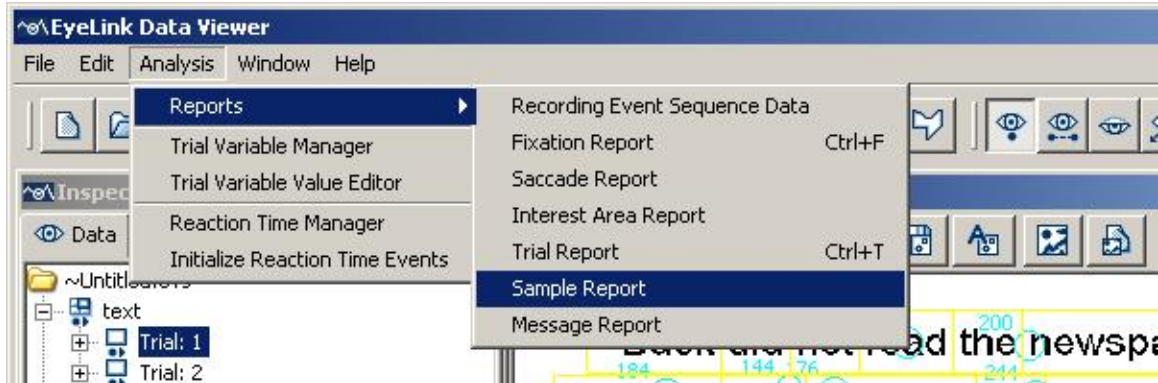


Figure 6-26. Creating a Sample Report.

2) In the following output report dialog, Select the variables in the list of available output variables (left panel) and press ">>" to enter the variables into the list of selected output variables.

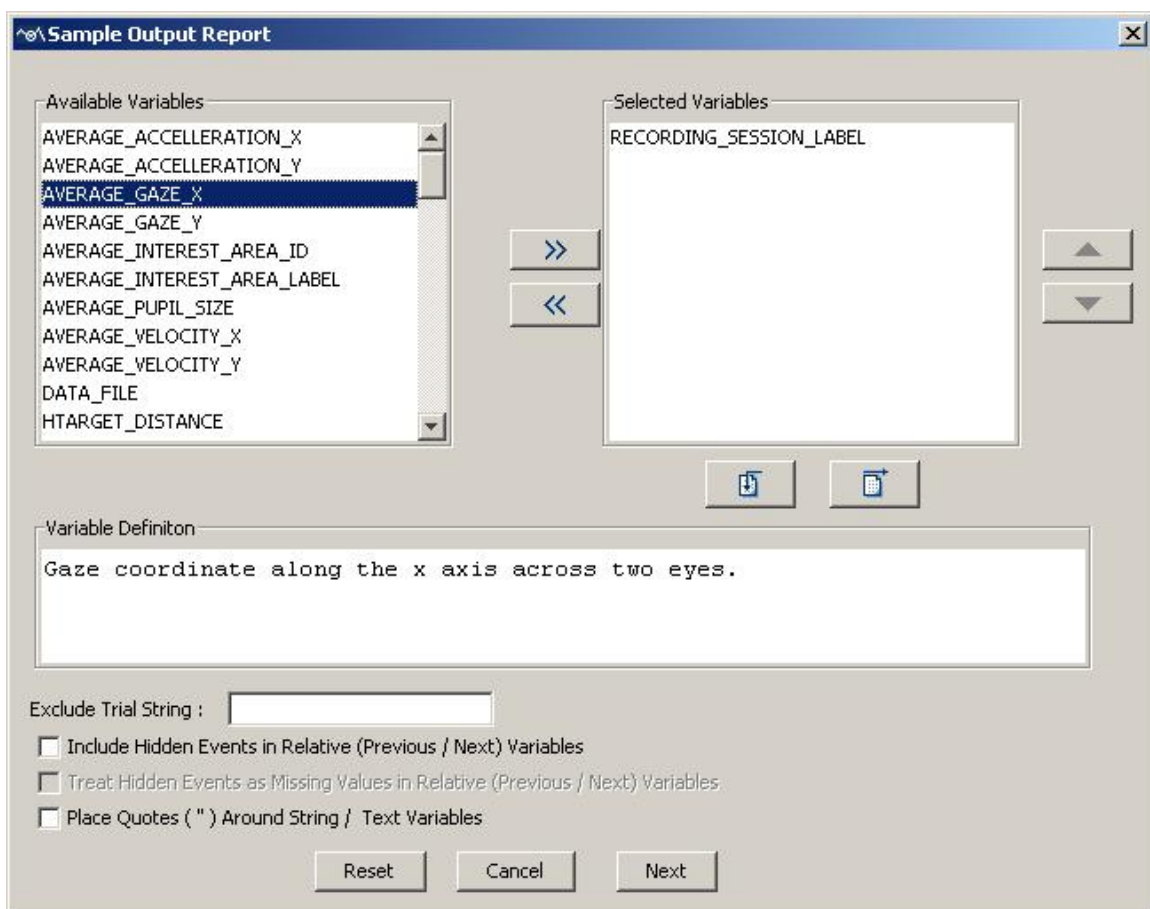




Figure 6-27. Variables in a Sample Report.

- To remove variables from the output list, simply select the variables and press <<.

- To change the order in which the variables are listed, simply select one variable and use the ▲ or ▼ button to move it to the desired position. Repeat this step until all variables are in the right position.
- Data Viewer supports exporting and importing of variable selection across viewing sessions. For example, the user may go through each of the output reports to configure the selection of variables in an intended order. Export the variable selection into a property file by clicking on the "Export Variable Selection" button . For all future viewing sessions, the variable selection in the reports can be automated by importing the previously saved property file (by clicking on the "Import Variable Selection" button ).

3) Once all the desired variables have been entered into the output list, press NEXT button.

4) In the Save dialog box, browse to the directory in which you want to save the output report. The default output directory can be configured through "General Preferences".

5) In the File name text box, enter the name for the document and choose the appropriate file extension. Note, if the output report file extension set to "All Files" and the file name does not contain an extension, the file name is appended with a ".txt" by default.

6) Press the "Save" button.

Note: Depending on the size of the EDF file(s), generating this report may take a long time.

6.9.3 Variables in the Sample Output Report

Currently, the sample output report provides output for the following variables:

Field	Contents
RECORDING_SESSION_LABEL	Label of the data file
AVERAGE_ACCELERATION_X	Acceleration value (in degree/second/second) along the x axis across two eyes ¹ .
AVERAGE_ACCELERATION_Y	Acceleration value (in degree/second/second) along the y axis across two eyes ¹ .
AVERAGE_GAZE_X	Gaze coordinate along the x axis across two eyes ¹ .
AVERAGE_GAZE_Y	Gaze coordinate along the y axis across two eyes ¹ .
AVERAGE_INTEREST_AREAS	List of interest areas in which the current sample (average across two eyes) falls.
AVERAGE_INTEREST_AREA_DA	Reports the data (interest area type,

TA	dynamic/static interest area, and coordinate data) of the interest area in which the current sample (average across two eyes) falls.
AVERAGE_INTEREST_AREA_DISTANCE	Distance (in degrees of visual angle) between the current sample (average across two eyes) and the center of the interest area in which the sample falls.
AVERAGE_INTEREST_AREA_ID	Index of the interest area in which the current sample (average across two eyes ¹) falls.
AVERAGE_INTEREST_AREA_LABEL	Label of the interest area in which the current sample (average across two eyes ¹) falls.
AVERAGE_INTEREST_AREA_PIXEL_AREA	Pixel area of the interest area in which the current sample (average across two eyes) falls.
AVERAGE_PUPIL_SIZE	Pupil size of the current sample across two eyes.
AVERAGE_VELOCITY_X	Velocity value (in degree/second) along the x axis across two eyes ¹ .
AVERAGE_VELOCITY_Y	Velocity value (in degree/second) along the y axis across two eyes ¹ .
DATA_FILE	File name of the recording;
HTARGET_DISTANCE	Distance between the head target and eye camera (in millimeters) for an EyeLink Remote recording. Returns a missing value if the head target was missing or if the data was recorded with a non-Remote eye tracker.
HTARGET_X	X position of the head target in camera coordinate for an EyeLink Remote recording. Returns a missing value if the head target was missing or if the data was recorded with non-Remote eye tracker.
HTARGET_Y	Y position of the head target in camera coordinate for an EyeLink Remote recording. Returns a missing value if the head target was missing or if the data was recorded with non-Remote eye tracker.
HTARGET_FLAGS	Flags used to indicate the head target tracking status ('.....' if target tracking is ok; otherwise error code) for an EyeLink Remote recording. Returns a missing value if the data was recorded with non-Remote eye tracker.
IP_DURATION	Duration of the interest period in milliseconds.
IP_END_TIME	End time (in milliseconds since EyeLink tracker was activated) of the interest period.

IP_LABEL	Label of the current interest period selected in the viewing session.
IP_START_TIME	Start time (in milliseconds since EyeLink tracker was activated) of the interest period. TRIAL_LABEL Label of the trial
LEFT_ACCELERATION_X	Acceleration value (in degree/second/second) of the left eye along the x axis.
LEFT_ACCELERATION_Y	Acceleration value (in degree/second/second) of the left eye along the y axis.
LEFT_FIX_INDEX	Ordinal sequence of the fixation in the trial to which the left-eye sample belongs.
LEFT_GAZE_X	Left eye gaze coordinate along the x axis.
LEFT_GAZE_Y	Left eye gaze coordinate along the y axis.
LEFT_INTEREST_AREAS	List of interest areas in which the left-eye sample falls.
LEFT_INTEREST_AREA_DATA	Reports the data (interest area type, dynamic/static interest area, and coordinate data) of the interest area in which the left-eye sample falls.
LEFT_INTEREST_AREA_DISTANCE	Distance (in degrees of visual angle) between the current left-eye sample and the center of the interest area in which the sample falls.
LEFT_INTEREST_AREA_PIXEL_AREA	Pixel area of the interest area in which the left-eye sample falls.
LEFT_INTEREST_AREA_ID	Index of the interest area in which the left-eye sample falls.
LEFT_INTEREST_AREA_LABEL	Label of the interest area in which the left-eye sample falls.
LEFT_IN_BLINK	Whether the left eye is in a blink.
LEFT_IN_SACCADE	Whether the left eye is in a saccade.
LEFT_PUPIL_SIZE	Left eye pupil size of the current sample.
LEFT_SACCADE_INDEX	Ordinal sequence of the saccade in the trial to which the left-eye sample belongs.
LEFT_VELOCITY_X	Velocity value (in degree/second) of the left eye along the x axis.
LEFT_VELOCITY_Y	Velocity value (in degree/second) of the left eye along the y axis.
RESOLUTION_X	Horizontal angular resolution (in screen pixels per degree) for the sample.
RESOLUTION_Y	Vertical angular resolution (in screen pixels per degree) for the sample.
RIGHT_ACCELERATION_X	Acceleration value (in degree/second/second) of

	the right eye along the x axis.
RIGHT_ACCELERATION_Y	Acceleration value (in degree/second/second) of the right eye along the y axis.
RIGHT_FIX_INDEX	Ordinal sequence of the fixation in the trial to which the right-eye sample belongs.
RIGHT_GAZE_X	Right eye gaze coordinate along the x axis.
RIGHT_GAZE_Y	Right eye gaze coordinate along the y axis.
RIGHT_INTEREST_AREAS	List of interest areas in which the right-eye sample falls.
RIGHT_INTEREST_AREA_DATA	Reports the data (interest area type, dynamic/static interest area, and coordinate data) of the interest area in which the right-eye sample falls.
RIGHT_INTEREST_AREA_DISTANCE	Distance (in degrees of visual angle) between the current right-eye sample and the center of the interest area in which the sample falls.
RIGHT_INTEREST_AREA_ID	Index of the interest area in which the right-eye sample falls.
RIGHT_INTEREST_AREA_LABEL	Label of the interest area in which the right-eye sample falls.
RIGHT_INTEREST_AREA_PIXEL_AREA	Pixel area of the interest area in which the right-eye sample falls.
RIGHT_IN_BLINK	Whether the right eye is in a blink.
RIGHT_IN_SACCADE	Whether the right eye is in a saccade.
RIGHT_PUPIL_SIZE	Right eye pupil size of the current sample.
RIGHT_SACCADE_INDEX	Ordinal sequence of the saccade in the trial to which the right-eye sample belongs.
RIGHT_VELOCITY_X	Velocity value (in degree/second) of the right eye along the x axis.
RIGHT_VELOCITY_Y	Velocity value (in degree/second) of the right eye along the y axis.
SAMPLE_BUTTON	If the EyeLink button box is pressed, this records the ID of the EyeLink button number.
SAMPLE_INDEX	The index of the sample in the trial.
SAMPLE_MESSAGE	Message text printed out during the current sample. Multiple messages associated with the current sample are delimited by a semicolon (;)
TARGET_ACCELERATION_X	Acceleration value (in degree/second/second) of the target along the x axis.
TARGET_ACCELERATION_Y	Acceleration value (in degree/second/second) of the target along the y axis.
TARGET_VELOCITY_X	Velocity value (in degree/second) of the target

	along the x axis.
TARGET_VELOCITY_Y	Velocity value (in degree/second) of the target along the y axis.
TARGET_VISIBLE	Whether the target is visible during the current sample. To support the retrieval of target position from a sample report, the target position data must be written according to some format (see section "Protocol for EyeLink Data to Viewer Integration -> Target Position Commands").
TARGET_X	Target gaze position on x axis.
TARGET_Y	Target gaze position on y axis.
TIMESTAMP	The time stamp of the sample (in milliseconds since EyeLink tracker was activated).
TRIAL_INDEX	Sequential order of the trial in the recording.
TRIAL_LABEL	Label of the trial
* TRIAL_START_TIME	The start time (in milliseconds since EyeLink tracker was activated) of the trial.
VIDEO_FRAME_INDEX	Index of the video frame that was visible at the sample time. Note that all "VIDEO_" variables are applicable to SceneLink recordings or some Experiment Builder recordings only. For SceneLink recordings, please check out the "Data Loading" Preference Settings.
VIDEO_NAME	Name of the video frame that was visible at the sample time.

Note ¹: All "AVERAGE_*" measures will have a missing value if both eye samples are not available (if pupils are missing or recording is done with left-eye or right-eye only).

In addition, other user-defined trial variable labels will also be listed. Please note that the data output in the sample report depends on the setting of the interest period filter (see section 6.4.2 or Figure 6-18) – only those fixation and saccade events falling within the active interest period will be reported. If a valid reaction time definition is found in a trial, the TRIAL_START_TIME will be set as the start as the reaction time definition; otherwise this will be set as the start of trial recording (see section 6.5.3 on discussion of three time periods used in Data Viewer).

6.10 Message Output Report

Message output report provides a columnar output of messages written in each trial as well as the eye movement events to which the message is associated. As with the other

reports, message report only outputs messages that fall within the active interest period for the trial.

6.10.1 Obtaining a Message Output Report

To obtain an output report for messages,

1) From the application menus, choose:

Analysis -> Report -> Message Report

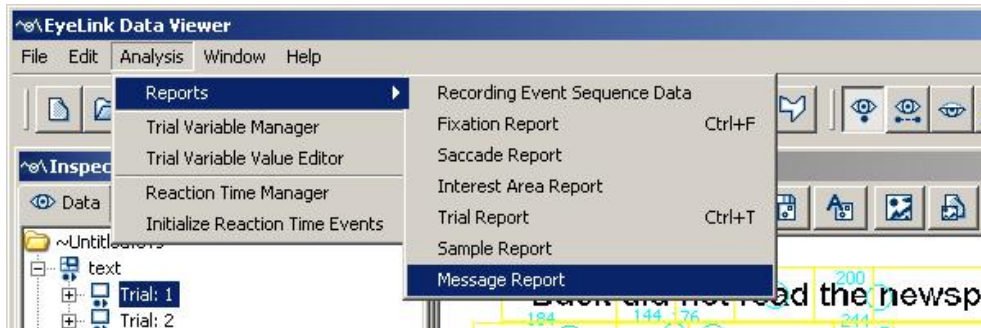


Figure 6-28. Creating a Message Report.

2) In the following output report dialog, Select the variables in the list of available output variables (left panel) and press ">>" to enter the variables into the list of selected output variables.

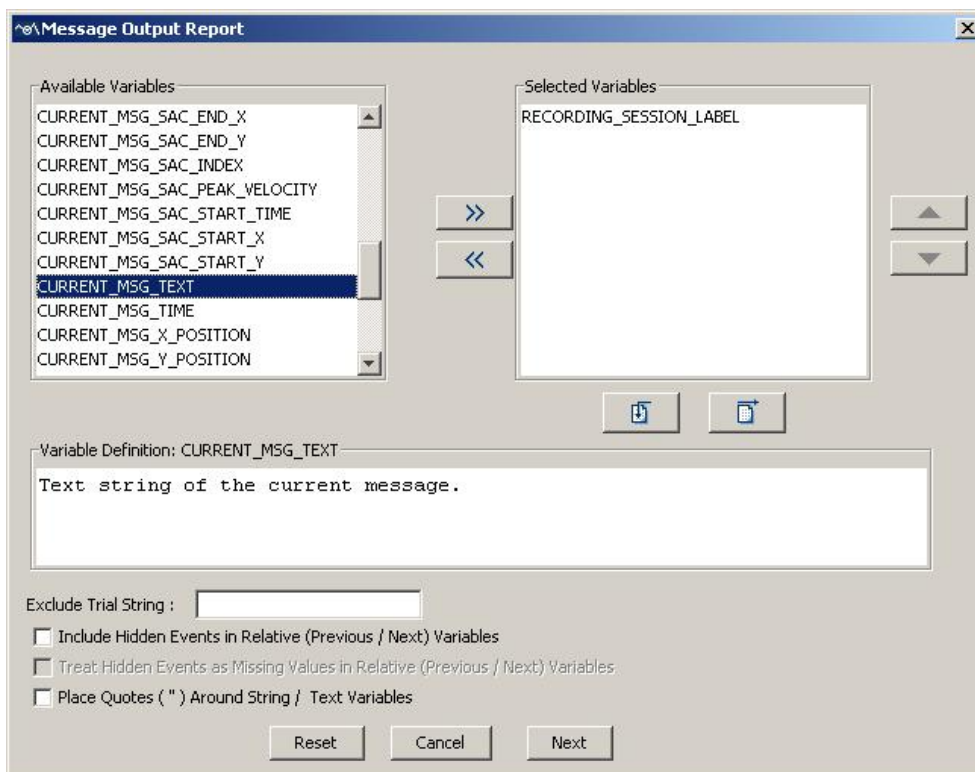




Figure 6-29. Variables in a Message Report.

- To remove variables from the output list, simply select the variables and press <<.
- To change the order in which the variables are listed, simply select one variable and use the ▲ or ▼ button to move it to the desired position. Repeat this step until all variables are in the right position.
- Data Viewer supports exporting and importing of variable selection across viewing sessions. For example, the user may go through each of the output reports to configure the selection of variables in an intended order. Export the variable selection into a property file by clicking on the "Export Variable Selection" button . For all future viewing sessions, the variable selection in the reports can be automated by importing the previously saved property file (by clicking on the "Import Variable Selection" button ).

3) Once all the desired variables have been entered into the output list, press NEXT button.

4) In the Save dialog box, browse to the directory in which you want to save the output report. The default output directory can be configured through "General Preferences".

5) In the File name text box, enter the name for the document and choose the appropriate file extension. Note, if the output report file extension set to "All Files" and the file name does not contain an extension, the file name is appended with a ".txt" by default.

6) Press the "Save" button.

Data Viewer also allows the user to configure the message output report. The user can exclude some trials by entering a string in the "Exclude Trial String" edit box. This string should be one message or part of a message within the scope of the trial recording that uniquely defines that subset of trials to be excluded. In addition, to enable correct loading of string/text variables in some statistics software, the Viewer also has an option of adding a pair of quotation marks for such variables.

6.10.2 Variables in the Message Output Report

Currently, the following variables are included in the message output report:

Field	Contents
RECORDING_SESSION_LABEL	Label of the data file
CURRENT_MSG_BLINK_DURATION	Duration of the blink to which the message is associated.

CURRENT_MSG_BLINK_END	End time (relative to the start of the interest period) of the blink to which the message is associated.
CURRENT_MSG_BLINK_INDEX	Index of the blink to which the message is associated.
CURRENT_MSG_BLINK_START	Start time (relative to the start of the interest period) of the blink during which the message was written.
CURRENT_MSG_FIX_DURATION	Duration of the fixation to which the message is associated.
CURRENT_MSG_FIX_END	End time (relative to the start of the interest period) of the fixation to which the message is associated.
CURRENT_MSG_FIX_INDEX	Index of the fixation to which the message is associated.
CURRENT_MSG_FIX_START	Start time (relative to the start of the interest period) of the fixation to which the message is associated.
CURRENT_MSG_FIX_X	X position of the fixation to which the message is associated.
CURRENT_MSG_FIX_Y	Y position of the fixation to which the message is associated.
CURRENT_MSG_INDEX	Index of the current message.
CURRENT_MSG_INTEREST_AREA	Interest areas in which the current eye position falls when the message is recorded.
CURRENT_MSG_INTEREST_AREA_ID	The ID of the interest area in which the current eye position falls when the message is recorded.
CURRENT_MSG_INTEREST_AREA_INDEX	The index of the interest area in which the current eye position falls when the message is recorded.
CURRENT_MSG_INTEREST_AREA_LABEL	Label for the interest area in which the current eye position falls when the message is recorded.
CURRENT_MSG_IS_RT_END	Whether the current message is the end event of a reaction-time definition.
CURRENT_MSG_IS_RT_START	Whether the current message is the start event of a reaction-time definition.
CURRENT_MSG_LABEL	Label of the current message event.
CURRENT_MSG_SAC_AMPLITUDE	Amplitude (in degrees of visual angle) of the saccade to which the message is associated.
CURRENT_MSG_SAC_AVG_VELOCITY	Average velocity (in degrees/second) of the saccade to which the message is associated.
CURRENT_MSG_SAC_DURATION	Duration of the saccade to which the message is associated.

N	associated.
CURRENT_MSG_SAC_END_TIME	End time (relative to the start of the interest period) of the saccade to which the message is associated.
CURRENT_MSG_SAC_END_X	X coordinate of the end point for the saccade to which the message is associated.
CURRENT_MSG_SAC_END_Y	Y coordinate of the end point for the saccade to which the message is associated.
CURRENT_MSG_SAC_INDEX	Index of the saccade to which the message is associated.
CURRENT_MSG_SAC_PEAK_VELOCITY	Peak velocity (in degrees/second) of the saccade to which the message is associated.
CURRENT_MSG_SAC_START_TIME	Start time (relative to the start of the interest period) of the saccade to which the message is associated.
CURRENT_MSG_SAC_START_X	X coordinate of the start point for the saccade to which the message is associated.
CURRENT_MSG_SAC_START_Y	Y coordinate of the start point for the saccade to which the message is associated.
CURRENT_MSG_TEXT	Text string of the current message.
CURRENT_MSG_TIME	Trial time (relative to the start of the interest period) of the message.
CURRENT_MSG_X_POSITION	X position of the current message text. It may be a missing value if samples are not loaded in the viewing session.
CURRENT_MSG_Y_POSITION	Y position of the current message text. It may be a missing value if samples are not loaded in the viewing session.
DATA_FILE	File name of the recording;
EYE_USED	Which eye's data (LEFT or RIGHT) is used to create this report?
IP_DURATION	Duration of the interest period in milliseconds.
IP_END_TIME	End time (in milliseconds since EyeLink tracker was activated) of the interest period.
IP_LABEL	Label of the current interest period selected in the viewing session.
IP_START_TIME	Start time (in milliseconds since EyeLink tracker was activated) of the interest period.
TRIAL_INDEX	TRIAL_LABEL Label of the trial Sequential order of the trial in the recording.
TRIAL_LABEL	Label of the trial

* TRIAL_START_TIME	The start time (in milliseconds since EyeLink tracker was activated) of the trial.
--------------------	--

Note: Messages written in the EDF file will be associated with an event closest in time. In very rare cases, the message can be written to both saccade and fixation events if the message was written out after the last sample of a fixation (or saccade) but before the first sample of the following saccade (or fixation).

In addition, other user-defined trial variable labels will also be listed. Please note that the data output in the message report depends on the setting of the interest period filter (see section 6.4.2 or Figure 6-18) – only those events falling within the active interest period will be reported. If a valid reaction time definition is found in a trial, the TRIAL_START_TIME will be set as the start as the reaction time definition; otherwise this will be set as the start of trial recording (see section 6.5.3 on discussion of three time periods used in Data Viewer).

6.11 Recording Event Sequence Data

Recording Event Sequence Data is a trimmed version of the EyeLink ASC file. For each trial, it outputs the selected events only and excludes the hidden events.

6.11.1 Obtaining a Recording Event Sequence Data

To obtain a recording event sequence data report:

- 1) From the menus, choose:
Analysis → Report → Recording Event Sequence Data.
- 2) In the Export dialog box, browse to the directory in which you want to save the data.
- 3) In the File name text box, type in the name for the data.
- 4) Press Export button.

6.11.2 Formats of the Recording Event Sequence Data

The following output is a trimmed version of a sample recording event sequence report.

```
MSG 5520005 DISPLAY_COORDS 0 0 1023 767
MSG 5520005 FRAMERATE 85.05 Hz.
MSG 5520009 TRIAL_VAR_LABELS PAGE
MSG 5684435 TRIALID PAGE1
START 5686598
BUTTON 5686573 5 true
MSG 5686595 DRIFTCORRECT R RIGHT at 320,40 OFFSET 0.09 deg. 4.1,1.4 pix.
```

```

MSG      5686597 RECCFG P 500 2 1
MSG      5686597 GAZE_COORDS 0.00 0.00 1023.00 767.00
EFIX     R 5686602 5686852 252 514.2 388.9 49.0   50.95 45.0   false [72: dogs, ]
BUTTON 5686716 5 false
MSG      5686723 DISPLAY ON
MSG      5686726 SYNCTIME 3
ESACC    R 5686854 5686900 48           514.1 385.7 342.1 84.5   7.58 229.0 50.55
         44.45
EFIX     R 5686902 5686980 80 340.2 80.6   47.0   50.2   43.9   false [5: the ]
ESACC    R 5686982 5687012 32 339.6 84.4   192.6 46.9   3.05 162.0 50.25 43.85
EFIX     R 5687014 5687176 164 195.0 44.4   46.0   50.3   43.8   false [3: not ]
ESACC    R 5687178 5687192 16 189.7 41.5   155.1 38.1   0.69 62.0 50.35 43.8
EFIX     R 5687194 5687364 172 153.8 45.2   44.0   50.4   43.8   false [2: did ]
ESACC    R 5687366 5687384 20 157.9 47.8   232.8 51.9   1.49 116.0 50.3   43.8
EFIX     R 5687386 5687588 204 233.9 54.8   45.0   50.25 43.8   false [3: not ]
ESACC    R 5718822 5718836 16 621.3 723.4   677.1 712.5 1.09 96.0 52.35 46.7
EFIX     R 5718838 5718968 132 684.3 710.4 46.0   52.45 46.7   false [142: sides.]
ESACC    R 5718970 5719032 64           685.0 706.9 347.0 106.4 14.87 371.0 51.3
         45.05
EFIX     R 5719034 5719336 304 356.5 131.3 43.0   50.25 43.5   false [16: not ]
BUTTON 5719243 5 true
MSG      5719244 ENDBUTTON 5
END 5719345
MSG      5719360 TRIAL OK

```

In this output file, the display coordinates and frame rate were reported first, followed by information from individual trials. This includes the selected fixations, saccades, blinks, messages, and button events and excludes any hidden events. For a Data Viewing session that loads in several recording sessions, only the first recording session will be exported in the recording event sequence output file.

6.11.2.1 Fixation Events

Fixation events are recorded in the format of:

- EFIX <eye> <stime> <etime> <dur> <axp> <ayp> <aps> <xr> <yr> <ma>
<ia>

This reports the eye used in the recording, the time of the first and last sample in the fixation, and computes the duration of the fixation in milliseconds. The average X and Y eye position (the type of position data is determined when the event was generated) and the average pupil size (area or diameter) are reported. The horizontal and vertical eye-position angular resolutions (in pixels per visual degree) are reported next. In addition,

the output contains information on whether the fixation has been manually adjusted <ma> and the ID and name of the interest area in which the current fixation falls <ia>.

6.11.2.2 Saccade Events

Saccade events are recorded in the format of:

- ESACC <eye> <stime> <etime> <dur> <sxp> <syp> <exp> <eyp> <ampl> <pv> <xr> <yr>

This reports the time of the first and last sample in the saccade, and computes its duration in milliseconds. The X and Y eye position at the start and end of the saccade (<sxp>, <syp>, <exp>, <eyp>) are listed. The total visual angle covered in the saccade is reported by <ampl>, which can be divided by (<dur>/1000) to obtain the average velocity. Peak velocity is given by <pv>. The horizontal and vertical eye-position angular resolutions (in pixels per visual degree) are given as well.

6.11.2.3 Blink Events

Blink events are recorded in the format of:

- EBLINK <eye> <stime> <etime> <dur>

This reports the eye used in the recording, the time of the first and last sample in the blink, and computes the duration of the blink in milliseconds.

6.11.2.4 Message Events

Messages are recorded in the format of:

- MSG <time> <message>

A message line contains the text of a time stamped message. The <message> text fills the entire line after the timestamp <time> and any blank space following it.

6.11.2.5 Button Events

Button events are recorded in the format of:

- BUTTON <time> <button #> <button_pressed>

Button lines report a change in state of tracker buttons. The <button #> reports which button has changed state. The <button_pressed> value will be true if the button has been pressed, false if it has been released.

7 Protocol for EyeLink Data to Viewer Integration

The EyeLink Data Viewer can interpret a set of message commands from an EyeLink data file that allow the viewer to automate some of the viewer configuration for a given data file or trial. Examples of these commands include defining the image to overlay and specifying trial variables. These messages can be written by the display experiment to the EDF file using the API function that allows the writing of a custom message to the EDF file.

Messages that make up the EyeLink data file protocol for the viewer are not added to the Message list that is created for viewing within the viewer tool. Message commands are therefore invisible to the viewing windows, and are only interpreted by the Viewer during application loading.

Important: Do not use space characters in any component of a viewer message command; the space character is used to tokenize and interpret all viewer messages. Instead of using a space character within a message token, use the _ character.

7.1 *Defining the Start and End of a Trial*

The viewer uses two special message commands to define what should be considered the start and end of a trial for the purpose of parsing the EyeLink data files.

7.1.1 Trial Start Message

Identifier (Default): TRIALID

Description:

This message defines the start of a trial for the viewer. This is different than the start of recording message START that is logged when the trial recording begins. The viewer will not parse any messages, events, or samples, that exist in the data file prior to this message. The command identifier can be changed in the data loading preference settings.

Format:

TRIALID <Trial ID value list>

Example:

MSG 3362269 TRIALID PIX1 images\sacrmeto.jpg

Note: If the trial start message is not found, or if no trial start message is defined for the viewing session, the START message is used to indicate the start of a trial for the Viewer.

7.1.2 Trial End Message

Identifier (Default): TRIAL_RESULT

Description:

This message defines the end of a trial for the viewer. This is different than the end of recording message END that is logged when the trial recording ends. The viewer will not parse any messages, events, or samples that exist in the data file after this message. The command identifier can be changed in the data loading preference settings.

Format:

TRIAL_RESULT <possible trial result values>

Example:

```
MSG 3383842 TRIAL_RESULT 0
```

Note: If the trial end message is not found, or if no trial end message is defined for the viewing session, the END message is used to indicate the end of a trial for the Viewer.

7.2 Pre-Trial Message Commands

The following message commands, if defined, MUST be sent to the EyeLink data file prior to the Trial Start message defined above (i.e., outside the scope of a trial as defined for the viewer). If they are sent during trial recording, these messages will not be interpreted.

7.2.1 Trial Variable Labels

Identifier (Default): TRIAL_VAR_LABELS

Description:

This message allows the definition of the labels to be used for trial variables within the viewer. The command identifier can be changed in the data loading preference settings.

Format:

TRIAL_VAR_LABELS <trial variable list>

Example:

```
MSG 3363553 TRIAL_VAR_LABELS TRIAL_VAR_1 TRIAL_VAR_2 TRIAL_VAR_3
```

7.2.2 Display Coordinates

Identifier: DISPLAY_COORDS or GAZE_COORDS

Description:

This message specifies the display coordinates to be used within the viewer. This message **MUST** be present for the viewer to function correctly.

Format:

```
DISPLAY_COORDS <left> <top> <right> <bottom>
```

Left, top, right, and bottom refer to the x-y coordinates of the top-left and bottom-right corners of display.

Example:

```
MSG 3325521 DISPLAY_COORDS 0 0 1279 1023
```

7.2.3 Trial Grouping

Identifier: V_TRIAL_GROUPING

Description:

Trial grouping message is used to group those trials in the same experimental condition together so that they can be manipulated (e.g., selecting, hiding, deleting, applying interest area templates) collectively. This message must be sent after TRIAL_VAR_LABELS message but before first trial start message in EDF file.

Format:

```
V_TRIAL_GROUPING <trial variable list>
```

The trial grouping variable(s) must be a member of the variables listed in the TRIAL_VAR_LABELS; otherwise, this variable is simply ignored.

Example:

```
MSG 3363553 V_TRIAL_GROUPING TYPE DIRECTION
```

Note: Trial variables are space delimited. Trial variables that do not exist in variable list for trial are ignored.

7.3 Trial Message Commands

The following message commands **MUST** be specified within the scope of a trial, or the message commands will not be interpreted by the viewer.

IMPORTANT: All message commands within the trial scope must start with a !V directive, or the command will not be interpreted and will instead appear as a regular message in the visible message list of the viewer.

Tip: In almost all instances, each of the following commands will be repeated for every trial of an EyeLink data file.

7.3.1 Trial Variable Values

Identifier (Default): TRIAL_VAR_DATA

Description:

This message specifies the list of trial variable values for the trial. The list of trial variable values must be in the same order as was specified in the TRIAL_VAR_LABELS command for the viewer to properly match the variable label with the variable value. The command identifier can be changed in the data loading preference settings.

Note: If this command is the same as the Trial Start Command (TRIALID), the values will still be extracted from the Trial Start Line; however a !V is not needed in the message.

Format:

```
!V TRIAL_VAR_DATA <trial_variable_list>
```

Example:

```
MSG 3363553 !V TRIAL_VAR_DATA value1 value2 value3
```

7.3.2 Single Trial Variable Message Token

Identifier (Default): TRIAL_VAR

Description:

This command lets the user specify a trial variable and value for the given trial. Send one message for each pair of trial condition variable and its corresponding value. This means that there is really no need to use TRIAL_VAR_LABELS command when using this command. The default command identifier can be changed in the data loading preference settings. Please note that the eye tracker can handle about 20 messages every 10 milliseconds. So be careful not to send messages too quickly when you have many trial condition messages to send. Add one millisecond delay between message lines if this is the case.

Format:

```
!V TRIAL_VAR <trial_var_label> <trial_var_value>
```

<trial_var_label> is the label for the variable being set. If the current Data Viewer session does not have this variable yet, it is created. <trial_var_value> defines the value to set for the variable for the given trial.

Example:

```
MSG 9350338 !V TRIAL_VAR trial 1
MSG 9350339 !V TRIAL_VAR condition gap
MSG 9350340 !V TRIAL_VAR direction Left
MSG 9350341 !V TRIAL_VAR gap_duration 200
MSG 9350342 !V TRIAL_VAR t_x 212
MSG 9350343 !V TRIAL_VAR t_y 384
```

7.3.3 Image Commands

The viewer supports a set of commands that display a default image on the overlay mode of the trial viewer screen. All image commands use the IMGLOAD token, followed by a sub command.

7.3.3.1 Image Loading – Fill Full Screen

Identifier: FILL

Description:

This message specifies the image to be used as the background for the spatial overlay view of a trial within the viewer. The image is sized to fit the dimensions specified in the DISPLAY_COORDS command message. The image should be represented as a relative path. The viewer will look for the image in the following order:

- 1) In the default image directory specified in the general preference settings.
- 2) In the directory the EyeLink data file is loaded from.
- 3) In the directory the viewer application is running from.

Format:

```
!V IMGLOAD FILL <relative_image_path>
```

Example:

```
MSG 3388468 !V IMGLOAD FILL Sac_blur.jpg
```

7.3.3.2 Image Loading – Top Left

Identifier: TOP_LEFT

Description:

This message specifies an image to be used as a segment of the spatial overlay view of a trial within the viewer. The image size is not changed (unless width and height parameters are also specified) and will be positioned by using the x and y positions specified as the top left corner of the image. The image should be represented as a relative path. The viewer will look for the image in the following order:

- 1) In the default image directory specified in the general preference settings.
- 2) In the directory the EyeLink data file is loaded from.
- 3) In the directory the viewer application is running from.

Format:

```
!V IMGLOAD TOP_LEFT <relative_image_path> <x_position>  
    <y_position> [width] [height]
```

Note: width and height are optional parameters

Example:

```
MSG 2740540 !V IMGLOAD TOP_LEFT fixations.gif 200 200
```

7.3.3.3 Image Loading – Centered

Identifier: CENTER

Description:

This message specifies an image to be used as a segment of the spatial overlay view of a trial within the viewer. The image size is not changed (unless width and height parameters are specified) and will be positioned centering on the x-y coordinate specified. The image should be represented as a relative path. The viewer will look for the image in the following order:

- 1) In the default image directory specified in the general preference settings.
- 2) In the directory the EyeLink data file is loaded from.
- 3) In the directory the viewer application is running from.

Format:

```
!V  IMGLOAD  CENTER  <relative_image_path>  <x_position>  <y_position>  
      [width]  [height]
```

Example:

```
MSG 2740540 !V IMGLOAD CENTER  fixations.gif 200 200
```

7.3.4 Simple Drawing

Besides the image loading, the viewer also supports a set of commands that draws simple graphics, such as lines, framed box, filled box, as well as display clearing.

7.3.4.1 Clear Overlay View

Identifier: CLEAR

Description:

This command clears the overlay view to the color specified by the red, blue, green integer values specified in the command. Each RGB value should be an integer between 0 and 255.

Format:

```
!V  CLEAR  <red>  <blue>  <green>
```


Example:

```
MSG 2740540 !V CLEAR 255 255 255
```

This clears the display to white color.

7.3.4.2 Line drawing

Identifier: DRAWLINE

Description:

This command draws a line with the color specified by the red, blue, green integer values specified in the command. Each RGB value should be an integer between 0 and 255.

Format:

```
!V DRAWLINE <red> <blue> <green> <x_start> <y_start> <x_end>  
          <y_end>
```

Note: <x_start>: x-coordinate of the starting point; <y_start>: y-coordinate of the starting point; <x_end>: x-coordinate of the end position; <y_end>: y-coordinate of the ending position.

Example:

```
MSG 9441901 !V DRAWLINE 255 0 0 100 100 300 400
```

This draws a red line, connecting (100, 100) and (300, 400).

7.3.4.3 Drawing a Rectangle

Identifier: DRAWBOX

Description:

This command draws a framed rectangle with the color specified by the red, blue, green integer values specified in the command. Each RGB value should be an integer between 0 and 255.

Format:

```
!V DRAWBOX <red> <blue> <green> <top> <left> <right> <bottom>
```

Note: Left, top, right, and bottom refer to the x-y coordinates of the top-left and bottom-right corners of rectangle.

Example:

```
MSG 9441901 !V DRAWBOX 0 255 0 200 300 400 500
```

This draws a green rectangle (frame), with the top-left corner at (200, 300) and the bottom-right corner at (400, 500).

7.3.4.4 Drawing a Filled Rectangle

Identifier: FILLBOX

Description:

This command draws a filled rectangle with the color specified by the red, blue, green integer values specified in the command. Each RGB value should be an integer between 0 and 255.

Format:

```
!V FILLBOX <red> <blue> <green> <top> <left> <right> <bottom>
```

Note: Left, top, right, and bottom refer to the x-y coordinates of the top-left and bottom-right corners of rectangle.

Example:

```
MSG 9441901 !V FILLBOX 0 0 255 400 500 500 600
```

This draws a rectangle, with the top-left corner at (400, 500) and the bottom-right corner at (500, 600). The rectangle is filled with the blue color.

7.3.4.5 Drawing a Fixation Point

Identifier: FIXPOINT

Description:

This command draws a fixation point at a specified position.

Format:

```
!V FIXPOINT <target_red> <target_blue> <target_green> <erase_red>
    <erase_green> <erase_blue> <x> <y> <outer_diameter>
    <inner_diameter>
```

Note: The RGB values for the fixation point drawing color is specified by the <target_red> <target_blue> <target_green> whereas the color of the fixation center is specified by <erase_red> <erase_green> <erase_blue>. The <outer_diameter> and <inner_diameter> specified the diameters of the drawing circles.

Example:

```
MSG 9441901 !V FIXPOINT 0 0 0 255 255 255 511 383 18 4
```

This draws a fixation point at (511, 383), with a diameter of 18 pixels for the outer ring and 4 pixels for the inner ring. The fixation point is drawn in black and its center is erased with white.

7.3.4.6 Draw List File

Identifier: DRAW_LIST

Description:

When a large number of simple drawings have to be made, it is easier to first record all of the simple drawings in a draw list file. During the viewing session, the viewer reads in the drawing commands from the pre-recorded file and re-creates the display drawing. In a draw list file (.dlf), each line represents one simple drawing or an image loading command written in the same format as mentioned above except that "!V" directive is not necessary.

Format:

```
!V DRAW_LIST <relative path>
```

Example:

```
Example: MSG 262495 !V DRAW_LIST BER\s1106.dlf
```

The content of the file being:

```
CLEAR 255 255 255
IMGLOAD TOP_LEFT BER\color.bmp 772 644 33 33
IMGLOAD TOP_LEFT BER\shape.bmp 772 314 33 33
IMGLOAD TOP_LEFT BER\orient.bmp 222 424 33 33
IMGLOAD TOP_LEFT BER\color.bmp 552 94 33 33
```

This first clears the display and then loads four small bitmap images.

7.3.5 Interest Area Commands

The viewer supports a set of interest area commands that inform the viewer what interest areas to create by default for a given trial. Trials that have interest area commands specified in the data file have a custom interest area set created when the data is loaded containing the specified interest areas.

All interest area commands start with the IAREA token and are followed by a one word sub command.

7.3.5.1 Rectangular Interest Area

Identifier (Default): RECTANGLE

Description:

This message specifies the attributes of a rectangular interest area for the trial. Each trial can have a set of such rectangular interest areas.

Format:

```
!V IAREA RECTANGLE <id> <left> <top> <right> <bottom> [label  
    string]
```

Note: The label string parameter is optional.

Example:

```
MSG 2740540 !V IAREA RECTANGLE 1 547 410 585 449 square
```

7.3.5.2 Elliptical Interest Area

Identifier (Default): ELLIPSE

Description:

This message specifies the attributes of an elliptical interest area for the trial. Each trial can have a set of such elliptical interest areas.

Format:

```
!V IAREA ELLIPSE <id> <left> <top> <right> <bottom> [label  
    string]
```

Note: The label string parameter is optional. Left, top, right, and bottom refer to the x-y coordinates of the top-left and bottom-right corners of the bounding rectangle.

Example:

```
MSG 2740540 !V IAREA ELLIPSE 2 647 310 685 349 circle
```

7.3.5.3 Freehand Interest Area

Identifier (Default): FREEHAND

Description:

This message specifies the attributes of a freehand interest area for the trial. Each trial can have a set of such freehand interest areas.

Format:

```
!V IAREA FREEHAND <id> < $x_1, y_1$ > < $x_2, y_2$ > ... < $x_n, y_n$ > [label string]
```

Note: The label string parameter is optional. x_n, y_n refers to the x, y coordinates of a point. The x, y coordinates of each point are separated by a comma.

Example:

```
MSG 2740540 !V IAREA FREEHAND 3 512,284 612,384 512,484 412,384 diamond
```

7.3.5.4 Interest Area Set

Identifier (Default): FILE

Description:

When a large number of interest areas have to be specified (such as in reading research), it is easier to first write all the interest area information in a text file and later on the Viewer can locate and import the interest-area file. The `!V IAREA FILE` serves this purpose. In an interest-area file, each line represents one interest area, recorded in the format of “RECTANGLE id left top right bottom [label]” for a rectangular interest area, “ELLIPSE id left top right bottom [label]” for an elliptical interest area, and “FREEHAND id x_1, y_1 x_2, y_2 ... x_n, y_n [label]” for a freehand interest area.

Format:

```
!V IAREA FILE <relative_file_path>
```

Example:

```
MSG 2740540 !V IAREA FILE segments\example.ias
```

7.3.6 Target Position Commands

For experiments in which a moving resource is used, the user may need to keep track of the position of the resource. The Data Viewer allows the user to calculate the position, velocity, and acceleration data for each individual samples in the sample output report. With such messages, position, velocity, and acceleration traces can be displayed in the time plot view.

Identifier (Default): TARGET_POS

Description:

This message specifies the position and visibility of the targets at the specific message time. The command identifier can be changed in the data loading preference settings. A semicolon is required if you are supplying two target positions.

Format:

```
!V TARGET_POS <Target1 Key> <(target1 x, target1 y)> <target 1  
visibility> <target 1 step>; <Target2 Key> <(target2 x, target2  
y)> <target 2 visibility> <target 2 step>
```

Where <Target1 Key> and <Target2 Key>: the tokens to extract the position(s) of target(s);
<(target1 x, target1 y)> and <(target2 x, target2 y)>: the position(s) of the targets. The coordinates must be enclosed in a pair of brackets;
<target 1 visibility> and <target 2 visibility>: if 1, the targets are visible; if 0, the targets are hidden;
<target 1 step> and <target 2 step>: if 1, the position of the target will be in steps; if 0, the position of the target is interpolated across samples.

Example:

```
MSG 9071000 -15 !V TARGET_POS TARG1 (351, 278) 1 0; TARG2 (590, 134)
1 0
```

Note: The user can use any token word to indicate the target ID. The "key" property of the temporal graph preference settings is set to blank by default so that various token words can be used. If two targets are involved and if the order of the two targets is important, set that field to the token strings used in the target position command string.

7.4 Reaction Time Definitions

To obtain an accurate reaction time (RT) measure, the EyeLink Data Viewer relies on a set of reaction time definitions to parse the start and end events for RT calculation. The reaction time definition command messages inform the Viewer of the RT start message, the type of RT end events (fixation, saccade, message, or button), the parameters of the end event, as well as the conditions to which the RT definition is applied. Reaction time definition commands can be provided in two ways. The user can employ a pre-trial (condition-based) command messages to specify a RT definition for each of the experimental condition. All pre-trial reaction time commands start with the V_RT token and are followed by a one word subcommand, indicating the type of RT end event (BUTTON, MESSAGE, FIXATION, or SACCAD). Alternatively, the user can supply a custom reaction time definition specifically for each trial. All custom trial-specific reaction time commands start with the !V V_CRT token and are followed by a one word subcommand.

Important: When an EyeLink Data File is loaded into the Viewer, any existing Reaction Time Definitions for the Data Viewing Session are removed and reloaded from the data file being imported. Therefore, it is important that each data file contain all the pre-trial (condition-based) command messages required for the full analysis, not just for the given data file. Any previously set custom, trial specific, reaction time definitions are not affected when a new data file is loaded.

7.4.1 Button Events

Identifier: BUTTON

Description:

This will treat a particular button press/release as the reaction time end event.

Pre-Trial Message Command Format:

```
V_RT  BUTTON  <rt_start_msg>  <button_id>  <state>  [CONDITION LABEL 1]
      [condition value 1]  [CONDITION LABEL 2]  [condition value 2] ...
```

Trial-specific Message Command Format:

```
!V  V_CRT  BUTTON  <rt_start_msg>  <button_id>  <state>
```

Note: Button_id refers to the intended button ID (from 1 to 8, use 0 for any button event). Button state could be 0 (button release) or 1 (button press). Condition labels and their values are not required for the trial-specific message command and optional for the pre-trial message command (if not specified, the RT definition is applied to all conditions). In case both a condition-based pre-trial message command and a custom trial specific message command are applicable for one trial, the trial-specific command will override the more general one.

Example:

```
MSG 329381 V_RT BUTTON SYNCTIME 5 1 TYPE normal
```

In this example, the end event for RT calculation is pressing button 5 serves. This RT definition is applicable to the condition in which the value of the variable TYPE is “normal”.

7.4.2 Fixation Events:

Identifier: FIXATION

Description:

This will end the trial by the appearance of a specific fixation. The user can specify the parameters of the fixation end event: the minimum fixation duration (in milliseconds) required, the *xy* coordinates for the region center, as well as the diameter of the fixation region.

Pre-Trial Message Command Format:

```
V_RT  FIXATION  <rt_start_msg>  <fixation_duration>  <x>  <y>
      <diameter>  [CONDITION LABEL 1]  [condition value 1]  [CONDITION
      LABEL 2]  [condition value 2] ...
```

Trial-specific Message Command Format:


```
!V V_CRT FIXATION <rt_start_msg> <fixation_duration> <x> <y>  
    <diameter>
```

Note: If the location information of the target fixation is not important, set the values of x , y , and diameter as -1. Condition labels and their values are not required for the trial-specific message command and optional for the pre-trial message command (if not specified, the RT definition is applied to all conditions).

Example:

```
MSG 329381 V_RT FIXATION SYNCTIME 500 -1 -1 -1
```

In this example, the RT end event is the occurrence of the first fixation with a minimum duration of 500 milliseconds whereas the location information of the fixation is not important.

7.4.3 Saccade Events:

Identifier: SACCADE

Description:

This will end the trial by the appearance of a specific saccade. The user can specify the parameters of the saccade end event: the minimum saccadic amplitude (in degrees of visual angle) required, the xy coordinates for the region center, as well as the diameter of the fixation region.

Pre-Trial Message Command Format:

```
V_RT SACCADE <rt_start_msg> <saccadic_amplitude> <x> <y> <diameter>  
    [CONDITION LABEL 1] [condition value 1] [CONDITION LABEL 2]  
    [condition value 2] ...
```

Trial-specific Message Command Format:

```
!V V_CRT SACCADE <rt_start_msg> <saccadic_amplitude> <x> <y>  
    <diameter>
```

Note: If the location information of the saccade is not important, set the values of x , y , and diameter as -1. Condition labels and their values are not required for the trial-specific message command and optional for the pre-trial message command (if not specified, the RT definition is applied to all conditions).

Example:

```
MSG 329381 V_RT SACCADE SYNCTIME 5.0 300 400 100 TYPE Normal
```

In this example, the RT end event is the occurrence of the first saccade with a minimum amplitude of 5.0 degrees of visual angle. The saccade must fall within 50 pixels from the region center (300, 400).

7.4.4 Message Events:

Identifier: MESSAGE

Description:

This will treat the first occurrence of a matching message as the end event for reaction time calculation.

Pre-Trial Message Command Format:

```
V_RT MESSAGE <rt_start_msg> <rt_end_msg> [CONDITION LABEL 1]  
[condition value 1] [CONDITION LABEL 2] [condition value 2] ...
```

Trial-specific Message Command Format:

```
!V V_CRT MESSAGE <rt_start_msg> <rt_end_msg>
```

Note: Condition labels and their values are not required for the trial-specific message command and optional for the pre-trial message command (if not specified, the RT definition is applied to all conditions).

Example:

```
MSG 329381 V_RT MESSAGE SYNCTIME END_RT
```

7.4.5 Reaction Time Definition Set

Identifier: FILE

Description:

When a large number of RT definitions have to be specified, it is easier to first record all RT definitions in a file and later on the viewer can read the RT definition file with the V_RT FILE command. In a reaction-time definition file, each line represents one RT definition, which is written in the same format as the pre-trial message command for each respective type of RT end event, except that in this case V_RT token is not necessary.

Pre-Trial Message Command Format:

V_RT FILE <relative path>

Example:

MSG 329381 V_RT FILE rt\default.rts

7.5 Changes in the Sample Experiments

The preceding sections list the special messages and commands that can be sent to the EDF file to enable the EyeLink Data Viewer to extract information in a more efficient way. The current section provides an overview of the changes made in the sample experiment templates. Please refer to the “Programming EyeLink Experiments in Windows Version 2.0” manual for the description of individual templates.

7.5.1 “Simple” Template

Two EyeLink Data Viewer commands have been introduced in the “Simple” template: “TRIAL_VAR_LABELS” for the definition of the labels to be used for trial condition variables within the viewer and “TRIAL_VAR_DATA” for the assignment of condition values for each trial. All these changes, listed in the following table, were made in the w32_simple_trials.c file of the template.

Data Viewer Functionality	Message Commands
Trial Variable Labels	eyemsg_printf("TRIAL_VAR_LABELS TRIAL_WORD");
Trial Variable Values	eyemsg_printf("!V TRIAL_VAR_DATA %s", trial_word[num-1]);

7.5.2 “Text” Template

Along with the “TRIAL_VAR_LABELS” and “TRIAL_VAR_DATA” commands, the “Text” template also uses “IMGLOAD FILL” to load a single image and “IAREA FILE” to import a file containing interest areas for the trial. All of these changes, listed in the following table, were made in the w32_text_trials.c file of the template.

Data Viewer Functionality	Message Commands
---------------------------	------------------

Trial Variable Labels	eyemsg_printf("TRIAL_VAR_LABELS PAGE");
Trial Variable Values	eyemsg_printf("!V TRIAL_VAR_DATA %d", num);
Images	eyemsg_printf("!V IMGLOAD FILL Images/%s", image_fn);
Interest Areas	eyemsg_printf("!V IAREA FILE Segments/%s", seg_fn);

7.5.3 “Picture” Template

One new feature of the “Picture” template is the demonstration of another approach to load images and to specify interest areas for the trial. The third trial of the template loads individual images (IMGLOAD TOP_LEFT) to create a composite image in the Viewer and reads individual rectangular interest areas directly (“IAREA RECTANGLE”). All these changes, listed in the following table, were made in the w32_picture_trials.c file of the template.

Data Viewer Functionality	Message Commands
Trial Variable Labels	eyemsg_printf("TRIAL_VAR_LABELS TYPE");
Trial Variable Values	eyemsg_printf("!V TRIAL_VAR_DATA %s", imgname[num-1]);
Images	
“Normal” image	eyemsg_printf("!V IMGLOAD FILL images/sacrmeto.jpg");
“Blurred” image	eyemsg_printf("!V IMGLOAD FILL images/sac_blur.jpg");
“Composite” image	eyemsg_printf("!V IMGLOAD TOP_LEFT %s %d %d %d %d", small_images[i], points[i].x, points[i].y, SCRWIDTH/2, SCRHEIGHT/2);
Interest Areas	
“Composite” image	eyemsg_printf("!V IAREA RECTANGLE %d %d %d %d %d %s", i+1, points[i].x, points[i].y, points[i].x + SCRWIDTH/2, points[i].y + SCRHEIGHT/2, small_images[i]);

7.5.4 “EyeData” Template

The following table lists all of the changes made to the w32_data_trials.c file of the “EyeData” template.

Data Viewer Functionality	Message Commands
Trial Variable Labels	eyemsg_printf("TRIAL_VAR_LABELS CONDITION");
Trial Variable Values	eyemsg_printf("!V TRIAL_VAR_DATA Playback");
Images	eyemsg_printf("!V IMGLOAD FILL images/grid.png");

Interest Areas	eyemsg_printf("!V IAREA FILE segments/grid.ias");
----------------	---

7.5.5 “GCWindow” Template

The following table lists all of the changes made to the w32_gcwindow_trials.c file of the “GCWindow” template.

Data Viewer Functionality	Message Commands
Trial Variable Labels	eyemsg_printf("TRIAL_VAR_LABELS TYPE CENTRAL PERIPHERAL");
Trial Variable Values	
Trial 1	eyemsg_printf("!V TRIAL_VAR_DATA TEXT TEXT MASK");
Trial 2	eyemsg_printf("!V TRIAL_VAR_DATA TEXT MASK TEXT");
Trial 3	eyemsg_printf("!V TRIAL_VAR_DATA IMAGE IMAGE MASK");
Trial 4	eyemsg_printf("!V TRIAL_VAR_DATA IMAGE MASK IMAGE");
Trial 5	eyemsg_printf("!V TRIAL_VAR_DATA IMAGE IMAGE BLURRED");
Images	
Trials 1 & 2	eyemsg_printf("!V IMGLOAD FILL images/text.png");
Trials 3, 4, & 5	eyemsg_printf("!V IMGLOAD FILL images/sacrmeto.jpg");
Interest Areas	
Trials 1 & 2	eyemsg_printf("!V IAREA FILE segments/text.ias");

7.5.6 “Control” Template

The following messages and commands have been added to the w32_control_trials.c file of the “Control” template.

Data Viewer Functionality	Message Commands
Trial Variable Labels	eyemsg_printf("TRIAL_VAR_LABELS CONDITION");

The following messages and commands have been added to the w32_control_trial.c file of the “Control” template.

Data Viewer Functionality	Message Commands
Trial Variable Values	eyemsg_printf("!V TRIAL_VAR_DATA GAZECTRL");
Images	eyemsg_printf("!V IMGLOAD FILL images/grid.png");
Interest Areas	eyemsg_printf("!V IAREA FILE segments/grid.ias");

7.5.7 “Dynamic” Template

One important feature introduced in the “Dynamic” template is the addition of custom reaction time definition (!V V_CRT”) for the saccade trials. Reaction time is defined as the interval between the onset of the target (“SYNCTIME” message) and the occurrence of the first saccade, with a minimum amplitude of 2.0°, that falls within 50 pixels from the saccade target (*goal_x*, *goal_y*). The following table lists all of the messages and commands that have been added in the w32_data_trials.c file of the “Dynamic” template.

Data Viewer Functionality	Message Commands
Trial Variable Labels	eyemsg_printf("TRIAL_VAR_LABELS TRIAL TYPE DIRECTION");
Trial Variable Values	eyemsg_printf("!V TRIAL_VAR_DATA %d %s %s", num, trial_labels[0][num-1], trial_labels[1][num-1]);
Reaction Time Definitions	
Trials 3-8	eyemsg_printf("!V V_CRT SACCAD E SYNCTIME 2.0 %d %d 50", goal_x, goal_y);

7.5.8 “Comm_simple” Template

Similar to the “Simple” template, “TRIAL_VAR_LABELS” and “TRIAL_VAR_DATA” commands were added to the w32_simple_trials.c file of the template (see the Table).

Data Viewer Functionality	Message Commands
Trial Variable Labels	eyemsg_printf("TRIAL_VAR_LABELS TRIAL_WORD");
Trial Variable Values	eyemsg_printf("!V TRIAL_VAR_DATA %s", trial_word[num-1]);



7.5.9 Other Templates

Since the “Broadcast” and “Comm_listener” templates do not produce EyeLink data recording, no change has been made in those templates to support Data Viewer analysis.

8 Preference Settings

All the elements of the EyeLink Data Viewer can be configured by the user from the Preference tab of the Inspector Window. These include everything from the default colors used, to the elements of how an EDF file is loaded. All the changes can be saved, allowing the changed preference settings to be loaded in subsequent viewing sessions.

Any preferences setting can be edited by clicking on the value cell and changing the value. For string and number values, be sure to press Enter after making the change to register your input with the application.

To make the changed preference settings persist over future sessions, the user can click the right mouse buttons and clicking on "Save Properties as Defaults". Data Viewer also supports exporting and importing of preference settings across viewing sessions. Export the preference settings into a property file can be done by clicking on the "Export properties" button . To use the saved preference settings for a given session, click on the "Import Properties" button .


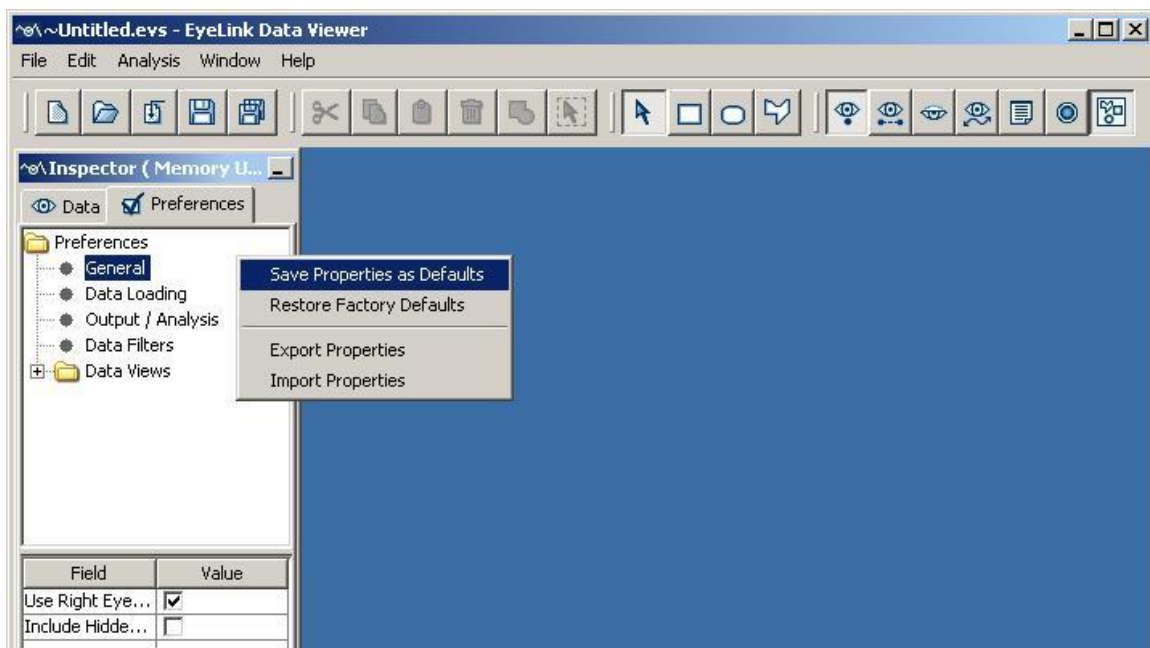


Figure 8-1. Saving properties as defaults.

8.1 General Preferences

The general preference setting covers the following elements:

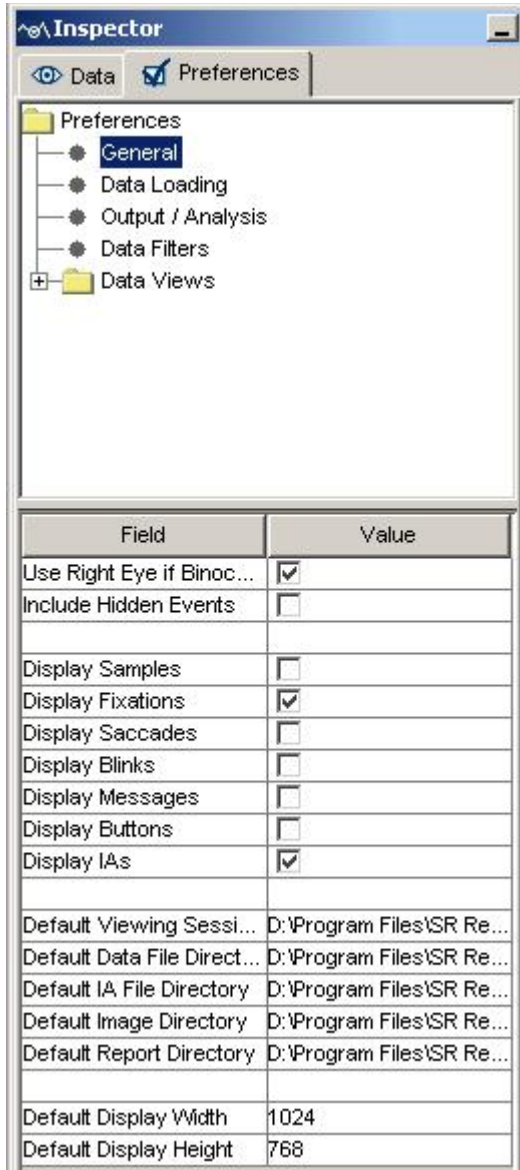

Use Right Eye if Binocular: For a binocular recording, which eye's data is to be displayed (right eye if checked; left eye if unchecked);

Include Hidden Events: Whether or not to include hidden events in Trial Views and Output Reports;


Display Samples, Display Fixations, Display Saccades, Display Blinks, Display Messages, Display Buttons, and Display IAs: Whether or not to display samples, fixations, saccades, blinks, messages, buttons, and interest areas when the data file is first loaded (by default, only fixations and interest areas are displayed);

Default Viewing Session File Directory, Default Data File Directory, Default IA File Directory, Default Image Directory, Default Video Directory, and Default Report Directory: Set default directories for accessing viewing session files, data file loading, interest area, images/video loading/saving, and data output. These settings can be changed by clicking on the current values and choosing a default directory from a Set window. Note, changes to these preference settings only take effect on files loaded after the change.

Default Display Width and Default Display Height: Set display width and height to enable the Data Viewer to run properly if DISPLAY_COORDS/GAZE_COORDS message is not found in the EDF file; otherwise, these preferences are ignored.

8.2 Data Loading Preferences

Data loading preference settings consist of the following parameters:

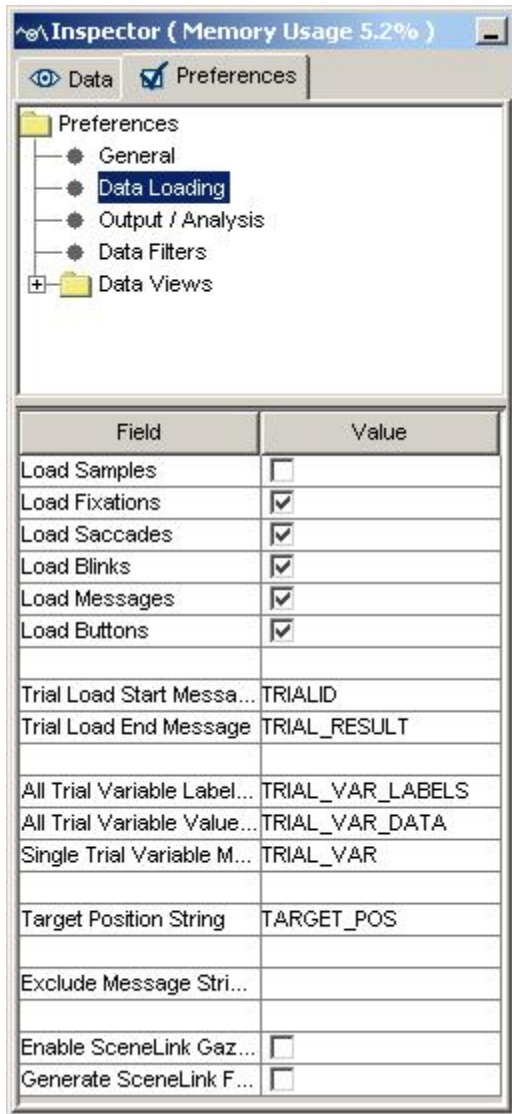

Load Samples, Load Fixations, Load Saccades, Load Blinks, Load Messages, Load Buttons: Allow the loading of samples, fixations, saccades, blinks, messages, and buttons from the EDF files. By default all of the events are loaded whereas samples are not. *Note:* Versions before 1.7.1 loads samples as well by default.

Trial Load Start Message and Trial Load End Message: Define the start and end messages of data loading. For each trial, all events/samples before the start message or after the end message will not be loaded. By default, the Viewer treats the “TRIALID” as the start message for data loading and “TRIAL_RESULT” as the end message. (Note: the **Trial Load Start Message** must be one message before START recording, otherwise the Viewer will internally re-set the loading position at the start of trial recording to keep trial data integrity.)

All Trial Variable Labels Message and All Trial Variable Values Message: The former allows the definition of labels to be used for trial variables within the viewer and the latter specifies the list of trial variable values for the trial. By default, the trial variable labels message is set as “TRIAL_VAR_LABELS” and the trial variable values message is set as “TRIAL_VAR_DATA”.

Single Trial Variable Message: This command lets the user specify a trial variable and value for the given trial. Send one message for each pair of trial condition variable

and its corresponding value. This means that there is really no need to use TRIAL_VAR_LABELS command when using TRIAL_VAR command. Default command token is "TRIAL_VAR".

Target Position String: String used to specify the position of the targets at the specific message time.

Exclude Message String: Any messages (within the scope of a trial as defined by the Trial Load Start Message and Trial Load End Message) containing the string will be ignored by the Viewer. Add a “;” to separate multiple message strings.

Collapse Identical Interest Areas with Different Time Stamps: If checked (the default setting), all interest areas with identical coordinates and label will be merged together (and thus treated as one single interest area). If unchecked, interest areas are organized based on the time stamps and therefore interest areas with identical coordinates and label but different time stamps are treated as separate items. In general, the former approach is typically used if the data analysis is done on the whole trial recording whereas the latter setting is used if the data analysis are done based on individual interest periods.

Enable SceneLink Gaze Mapping: If checked, allows the user to load EDF file with mapped gaze data to scene video coordinates.


Generate SceneLink Frame Message: If checked, allows the user to generate frame number messages when loading file. The frame number message will only be written if a lookup table has been created for synchronized playback of eye movement data over the scene video (see section 5.1 of the "EyeLink II Scene Camera User Manual").

Default Encoding: This option is used in case the interest area files are saved in a different encoding type.

Note: The user has to save the preferences and reload the data file before the changes in the above preference settings take effect.

8.3 Output / Analysis

The output / analysis preference settings consist of the following elements:

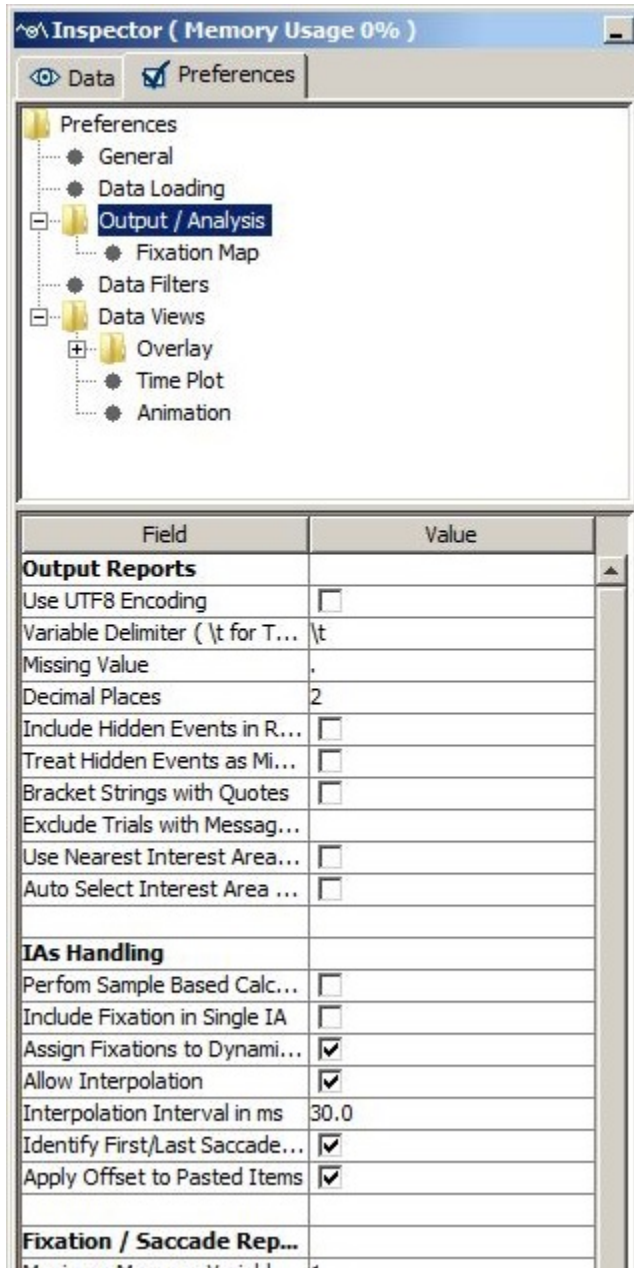

Output Reports

Use UTF8 Encoding: By default (unchecked), the saved output file will be a plain text file in ANSI encoding. If checked, the saved tab-delimited text file will be in UTF-8 encoding and the .XLS file will be in Unicode Text to better support non-ASCII characters in the output reports.

Variable Delimiter (\t for Tab):

Specifies the delimiter string for separating columns in the output file.

Missing Value: Replaces variables with missing value with a string specified here.

Decimal Places: Specifies the number of digits to appear after the decimal point.

Include Hidden Events in Relative Variables: Allows the inclusion of hidden events in the relative event variables [i.e., those variables beginning with PREVIOUS_FIX_, NEXT_FIX_, PREVIOUS_SAC_, or NEXT_SAC_].

Treat Hidden Events as Missing Values: If hidden events in the relative variables are included in the analysis, the user can further decide whether or not to treat these events as missing data.

Bracket Strings with Quotes: Adds a pair of quotation marks for string variables.

Exclude Trials with Message String: Excludes trials containing the specified message string. This string should be one message or part of a message within the scope of a trial that uniquely defines that subset of trials to be excluded.

Auto Select IAS for Trial: If true, trial interest area set association is done automatically by using "Image Name mapping" attribute of interest area set templates in the Data Viewer session (see section 5.8.8 "Automatic Trial Image-Interest Area Set Association" for details). The default is false.

Use Nearest Interest Area for Outlier Fixations: If the current fixation is not in any defined interest areas, assigns the nearest interest area as the current interest area.

IAs Handling:

Perform Sample Based Calculation: If selected, samples within fixations (if loaded) are used for interest area matching and report calculation. If disabled (default option), the average x, y position of the fixation is used. This option makes a difference for those data sets where fixations have a large spatial spread (e.g., smooth pursuit).

Include Fixation in Single IA: When selected, the fixation is assigned to one of the interest areas only - the interest area assigned gets the full length of the fixation when calculating dwell times. If this option is disabled (default setting), the fixation can be assigned to multiple interest areas, and effective duration (the total amount of time that fixation and instances of the same IA intersect) for each of the interest areas is used for the calculation of dwell times.

Assign fixations to dynamic IA ordered by relative duration: Some variables (e.g., "CURRENT_FIX_INTEREST_AREA_INDEX" in the Fixation Report) require a fixation to be assigned to a single interest area. If a fixation lands in multiple dynamic interest areas, the assignment by default will be based on the duration of fixation on respective interest areas - the interest area with the longest duration gets the fixation assignment. Unchecking this option will do the fixation assignment by interest area index.

Allow Interpolation: If selected, any dynamic interest area that has its interpolation property enabled will interpolate specially across its duration to the next instance's position and shape (if it's of the same shape type).

Interpolation Interval in ms: How often are dynamic interest areas interpolated?

Identify First/Last Saccade Based on First/Last Fixation: The end point of a saccade doesn't necessarily have the same coordinate as the following fixation event (because the eye is not absolutely still during a fixation; small eye movements such as drifts, tremors, microsaccades can occur during a fixation). As a result, a saccade and the following fixation may end into different interest areas. If the "Identify First/Last Saccade Based on First/Last Fixation" option is checked, Data Viewer first identifies the first and last fixation event in a trial and then find the preceding saccade of the fixation event and use this as the first/last saccade of the trial, despite occasionally the identified saccade may not end into the current interest areas. If this option is unchecked, the saccades are identified based on the actual end point and there might be a disconnect between the first/last saccade and the first/last fixation identified for the trial.

Apply Offset to Pasted Items: If checked, then pasted interest areas or interest area instances will be offset slightly down and to the right from the location of the original interest area or interest area instance that has been cut/copied to the clipboard.

Fixation/Saccade Report


Maximum Message Variables for Eye Event: Maximum number of messages associated with a given fixation or saccade that will be shown in the output report.

Eye Event Msg. Variable String Match Token: If left blank, all messages associated with an eye event will be reported; otherwise, only those messages containing the "token" will be reported.

Fixation Stream Analysis

Fixation Skip Count: Defines the number of fixations to skip when looking for the next interest area to use in the fixation sequence analysis variables of the Interest Area report. Default is 1. Cannot be less than 1.

8.4 Fixation Map

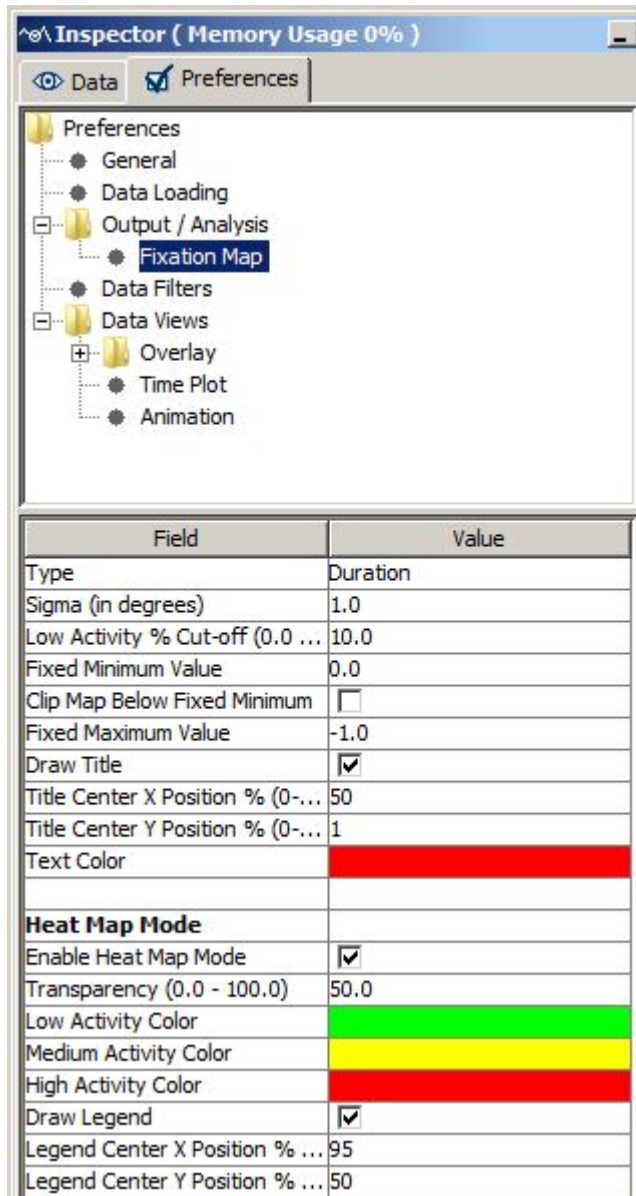

Type (1 = Duration, 2 = Count, 3 = Count Density, 4 = Duration Density):

The fixation map is created based on the duration of fixations (1), count of the fixation (2), probability of fixation across the display (3), or proportion of trial dwell time (4).

Sigma (in degrees): Sets the standard deviation of the Gaussian distribution for each fixation point when creating a fixation map.

Low Activity % Cut-off (0.0 - 100.0): Sets the minimum amount of activation required to have an area painted with an activation color - this is a certain percent of the actual maximum per-trial-average value for the map to keep the contour of the fixation map in case the actual maximum per-trial-average value is different from the "Fixed Maximum Value". If set to 0, the whole map is painted (with most of the areas in low-activity color).

Fixed Minimum Value: By default, the scale of the legend starts from 0. A non-zero value set here will be used as the minimum value of the scale. Any areas with an activation value below this "fixed minimum value" but above the "Low Activity % Cut-off (0.0- 100.0)" threshold, will be painted with the low activity color.

Clip Map below Fixed Minimum: If this enabled, this will not draw an areas with an activation value below the "Fixed Minimum Value" set above.

Fixed Maximum Value: If set to the default value of -1, the maximum value in the legend is based on the actual heat map maximum value. Any value > 0 will be used as the maximum legend scale. Any areas with an activation value above this "fixed maximum value" will be painted with the high activity color.

Draw Title: If checked, a title is displayed on the fixation map.

Title Center X Position % (0-100): The X position for the center of the title in percent of the display width.

Title Center Y Position % (0-100): The Y position for the center of the title in percent of the display height.

Text Color: Sets the color for the text of the heat map title and legend.

Enable Heat Map Mode: If checked, the fixation map will be drawn in a colored heat map; if unchecked, a grayscale version will be created.

Transparency (0.0 - 100.0): Sets the transparency level of the heat map. If 100, the heat map will be solid.

Low Activity Color: Sets the color used to draw low activity area.

Medium Activity Color: Sets the color used to draw medium activity area.

High Activity Color: Sets the color used to draw high activity area.


Draw Legend: If checked, a color scale is displayed on the fixation map.

Legend Center X Position % (0-100): The X position for the center of the legend in percent of the display width.

Legend Center Y Position % (0-100): The Y position for the center of the legend in percent of the display height.

8.5 Data Filter Preferences

The data filter preference settings consist of the following elements:

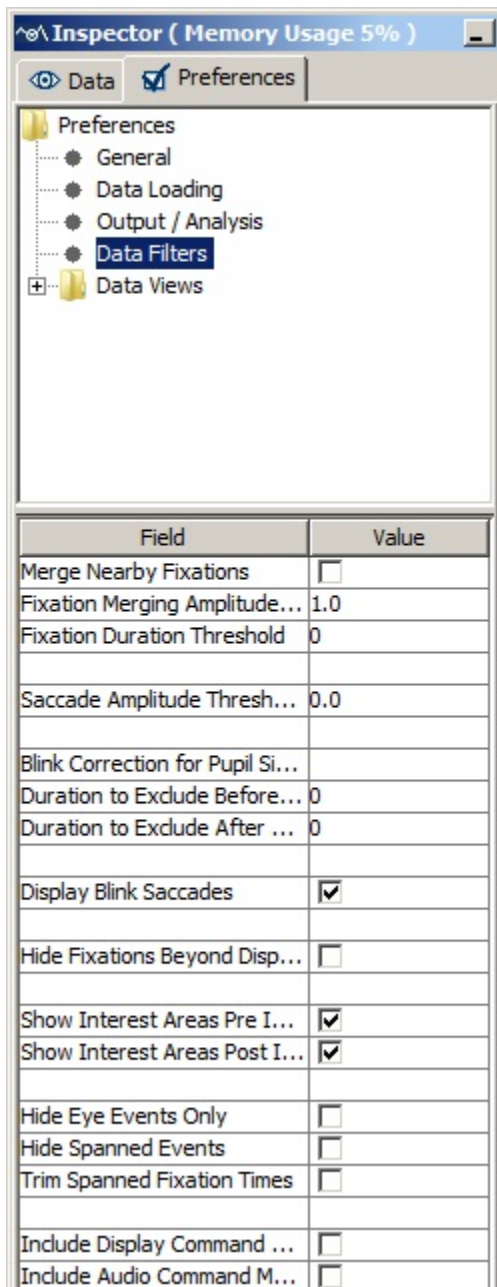

Merge Nearby Fixations: If checked, merge the neighboring brief fixations. A fixation to be merged must have a duration of less than that specified by Fixation Duration Threshold and be within 1° of the target fixation. These default thresholds are set in the “Amplitude Threshold” and “Fixation Duration Threshold” fields.

Fixation Merging Amplitude Threshold: Sets the amplitude threshold for fixation merging. The default value is 1.0°.

Fixation Duration Threshold: Sets the fixation duration threshold for fixation merging. A fixation will be merged if “Merge Nearby Fixations” is on and its duration is less than the specified duration threshold.

Saccade Amplitude Threshold: Sets the minimum size of a saccade to be displayed in the viewing session. All saccades will be displayed with an amplitude threshold set to 0.0.

Blink Correction for Pupil Size Calculation: Duration to Exclude Before Blinks (msec): Milliseconds of samples to exclude before the start of the blink events for pupil size calculation.

Duration to Exclude After Blinks (msec): Milliseconds of samples to exclude after the end of the blink events for pupil size calculation.

Display Blink Saccades: Shows flanking saccades of a blink.

Hide Fixations Beyond Display Bounds: Whether or not to hide the fixations falling beyond the display boundaries;

Show Interest Areas Pre Interest Period: If interest periods are defined, whether or not to display interest area timed before the start of a selected interest period. If no interest periods are defined, the interest areas are displayed anyway.

Show Interest Areas Post Interest Period: If interest periods are defined, whether or not to display interest area timed after the end of a selected interest period. If no interest periods are defined, the interest areas are displayed anyway.

Hide Eye Events Only: If checked, filters out eye events (fixations, saccades, and blinks) falling outside of the RT definition. If unchecked, filter out all events, including buttons events and messages.

Hide Spanned Events at Start of IP: Hide fixations that start before the beginning of the currently-selected interest period and end after the beginning of the currently-selected interest period. This will hide them from the various views of Data Viewer and will exclude them completely from Reports and Fixation Maps.

Hide Spanned Events at End of IP: Hide fixations that start before the end of the currently-selected interest period and end after the end of the currently-selected interest period. This will hide them from the various views of Data Viewer and will exclude them completely from Reports and Fixation Maps.

Trim Spanned Fixation Duration at Start of IP: Trim the duration of fixations that start before the beginning of the currently-selected interest period and end after the beginning of the currently-selected interest period. Trimming them will effectively exclude the portion of the spanned fixation that falls outside the interest period. This will influence the spanned fixations' durations in the various views of Data Viewer and in the output that goes into Reports and Fixation Maps.


Trim Spanned Fixation Duration at End of IP: Trim the duration of fixations that start before the end of the currently-selected interest period and end after the end of the currently-selected interest period. Trimming them will effectively exclude the portion of the spanned fixation that falls outside the interest period. This will influence the spanned fixations' durations in the various views of Data Viewer and in the output that goes into Reports and Fixation Maps.

Include Display Command Messages: Whether display command messages (i.e., command message starts with “!V” listed in sections 7.3.2 image loading and 7.3.3 simple graphics) should be visible.

Include Audio Command Messages: Whether audio command messages should be visible. The following built in audio messages are supported: "!V ARECSTART", "!V ARECSTOP", "!V VOICEKEY", "!V APLAYSTART", and "!V APLAYSTOP".

8.6 General Trial View Preferences

The parameters related to the general trial view preference settings are covered below:

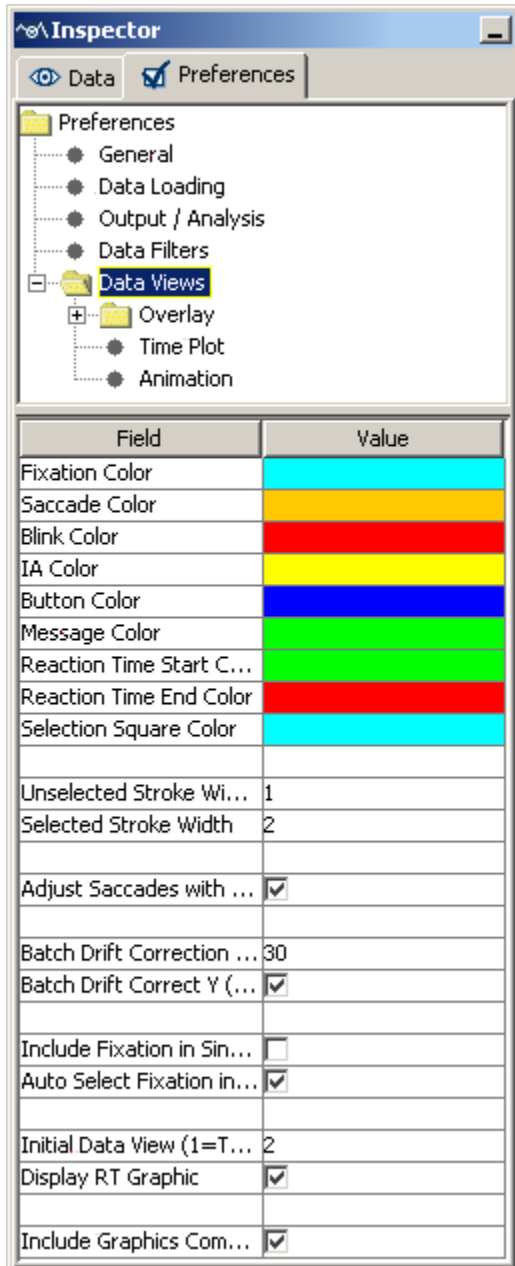

Fixation Color, Saccade Color, Blink Color, IA color, Button Color, Message Color, Reaction Time Start Color, Reaction Time End Color, and Selection Square Color: Set the colors to be used to show fixations, saccades, blinks, interest areas, button events, messages, the start/end of reaction time calculation, and the event selection square. These settings can be changed by clicking the corresponding cell and choosing a color from the palette dialog or entering the value in RGB or HSB scales.

Unselected Stroke Width and Selected Stroke Width: Pixel width for the selected and unselected events;

Adjust Saccades with Fixations: Whether or not to adjust saccade parameters when the fixations are shifted or merged;

Batch Drift Correction Error Threshold: The maximum tolerable deviation in the Y dimension (or the X dimension if the next field is unchecked) for a fixation from the mean of the group of fixations to be drift-corrected in a batch. If any fixations within the group exceeds this threshold, batch drift correction cannot be performed.

Batch Drift Correction Y (X=false): If checked, chooses the Y dimension for batch drift correction. If unchecked, X dimension is checked;

Auto Select Fixation in Selected IA: When an interest area is selected, whether or not to select fixations inside the IA as well;


Initial Data View (1 = TP; 2 = OL; 3 = AV): Selects the mode (1= temporal graph mode; 2 = spatial overlay mode; 3 = animation playback mode) of the trial view window when the data is first loaded;

Display RT Graphics: If reaction time definitions are initialized, whether or not to show the “RT” symbol in the spatial overlay view and to draw the Reaction Time Start/End lines in the temporal graph view of the trial data.

Include Graphics Commands Sent Prior to IP: If no graphics commands are found within the current interest period, uses those sent prior to the current interest period.

8.7 Spatial Overlay View Preferences

The spatial overlay view preference settings cover the following parameters:

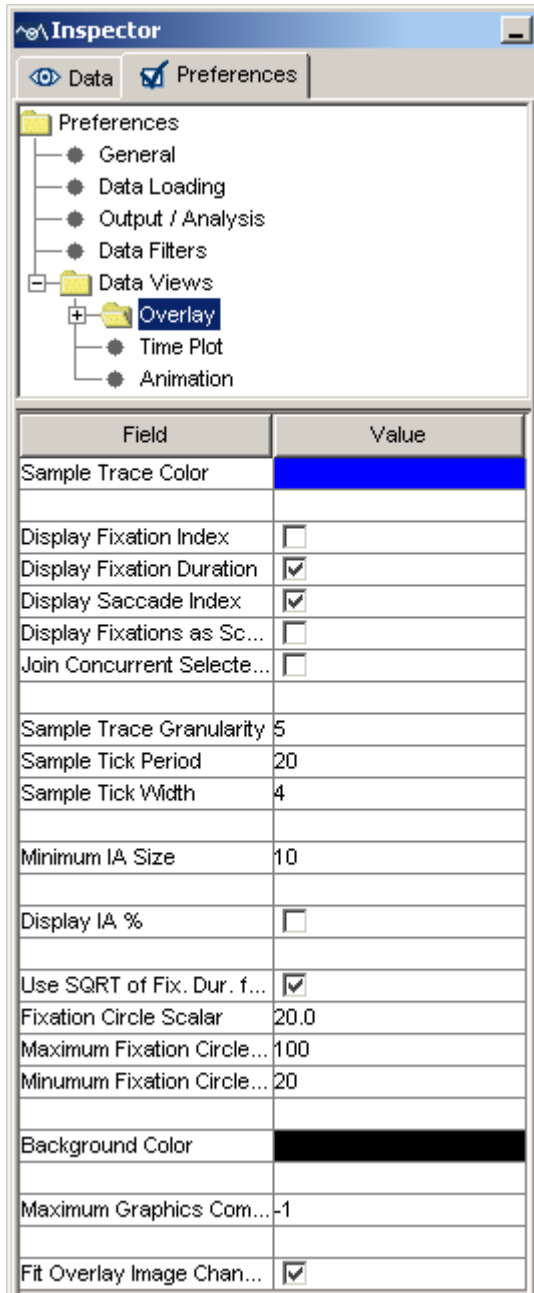

Sample Trace Color: Sets the color to be used to show the sample trace;

Display Fixation Index, Display Fixation Duration, and Display Saccade Index: Whether or not to show the fixation index/fixation duration/saccade index in the spatial overlay view;

Display Fixations as Scan Path: If checked, two temporally contiguous fixations are connected by an arrow, pointing towards the more recent fixation;

Joint Concurrent Selected Fixations: For the selected fixations, connects temporally contiguous fixations;

Display Text Font Size: Size of the font used in the Overlay view (e.g., size of the font for fixation durations/indices).

Sample Trace Granularity: If the sample visibility is toggled on, shows every n th sample (by default, $n = 5$);

Sample Tick Period: Draws a circular tick for every n th sample (by default, $n = 20$). Widely spread ticks indicate high saccade velocity;

Sample Tick Width: Diameter (in pixels) of the sample ticks;

Minimum IA Size: Minimum width of an interest area;

Display IA%: If selected, shows percentage of trial dwell time spent on the selected interest area

Use SQRT of Fixation Duration for Circle Size: Uses the square root of fixation duration to calculate the width of the fixation circle if checked.

Fixation Circle Scalar, Maximum Fixation Circle Width, and Minimum Fixation Circle Width: If the above **Use SQRT of Fixation Duration for Circle Size** option is not used, use **Fixation Circle Scalar** to define a scalar for fixation circle width calculation. The maximum and minimum circle widths are set with the **Maximum Fixation Circle Width/Minimum Fixation Circle Width** elements;


Background Color: Background color for the spatial overlay view (used only if the trial bitmap is not loaded).

Maximum Graphics Commands to Process: Sets the maximum number of graphics commands to process in each trial. This command is useful for those trials in which there are many viewer graphics commands while the user may just need to see a few of them. If it is set to -1, the viewer will process all graphics commands. For all other entries of integer numbers, the viewer will process the number of graphics commands specified when creating background image for overlay view.

Fit Overlay Image Change to Display: If True, when the background image for a trial is manually changed, the image is scaled to fit the display dimensions for the trial. If false, the image size is not altered and the image is centered on the trial overlay view.

8.8 Image Segmentation Preferences

The parameters related to the auto image segmentation preference settings are covered below:

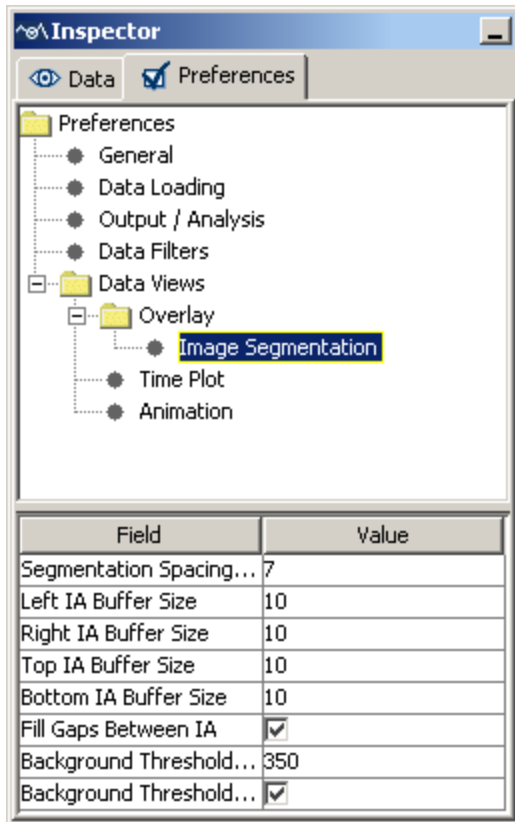

Segmentation Spacing Threshold: Number of consecutive pixels below threshold before segment end is identified.

Left IA Buffer Size, Right IA Buffer Size, Top IA Buffer Size, Bottom IA Buffer Size: Number of pixels added to the left, top, right, and bottom of the interest area. Note that the **Top IA Buffer Size, Bottom IA Buffer Size** fields will not have an effect if **Fill Gaps Between IA** is checked.

Fill Gaps Between IA: If checked, gaps between consecutive Interest Areas will be filled by expanding the size of each Interest Area.

Background Threshold (0-765): Combined RGB (R+G+B) to be used as the threshold limit for segmentation.


Background Threshold Less Than: If checked, pixels with a RGB value less than the threshold color are considered 'background' pixels. This setting is good for displays with a lighter background. If this field is unchecked, pixels with an RGB value greater than the threshold are considered background pixels (good for displays with a darker background).

Segmentation Direction: If "Left to Right" is selected then the interest areas will be created such that the interest areas furthest to the left will have the lowest index/id and the index/id values will increase in a left-to-right manner. If "Right to Left" is selected then the interest areas will be created such that the interest areas furthest to the right will have the lowest index/id and the index/id values will increase in a right-to-left manner.

Note: Auto segmentation may not work for a display with vertical stripes along its border(s) that have the 'foreground' color. If such stripes are not part of the original image, try to reload the background image or to change the background color for the spatial overlay view to match with the background color of the original image.

8.9 Temporal Graph Preferences

The parameters related to the temporal graph view preference settings are covered below:

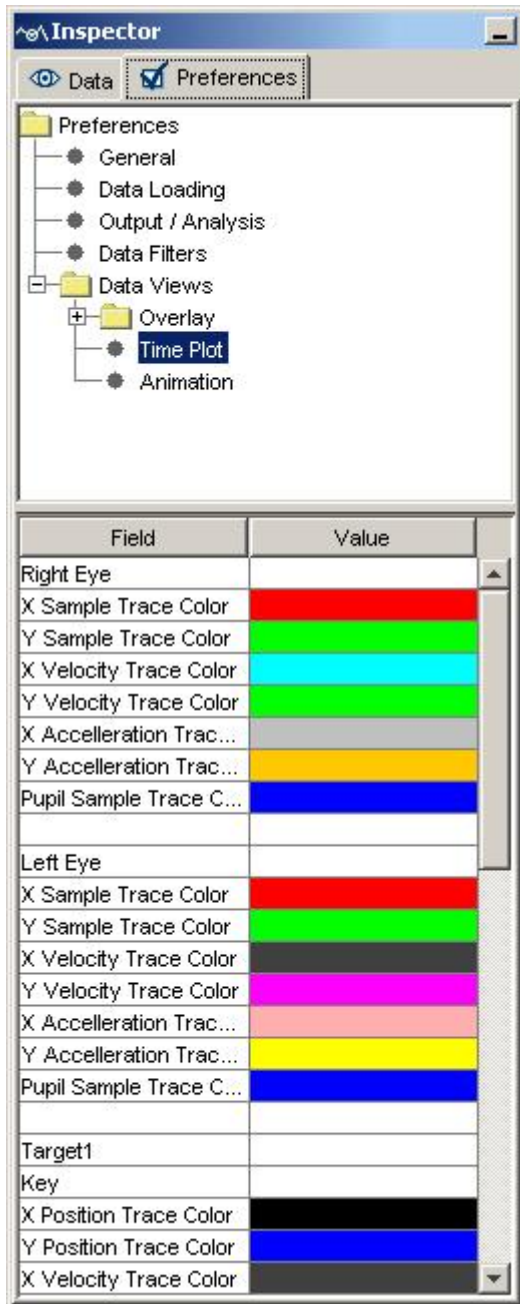

Left Eye/Right Eye

X Sample Trace Color, Y Sample Trace Color, X Velocity Trace Color, Y Velocity Trace Color, X Acceleration Trace Color, and Pupil Sample Trace Color: Set the colors to be used to show the horizontal and vertical position, velocity, and acceleration traces of the eye samples. The pupil sample traces of both eyes are also displayed. The default settings can be changed by clicking the corresponding cell and choosing a color from the palette or entering the value in RGB or HSB scales;

Target 1/Target 2

Key: Token word to indicate the target ID. This field is set to blank by default so that various token words can be used. If two targets are involved and if the order of the two targets is important, set this field to the token string used in the EDF file.

X Position Trace Color, Y Position Trace Color, X Velocity Trace Color, Y Velocity Trace Color, X Acceleration Trace Color, and Pupil Sample Trace Color: Set the colors to be used to show the horizontal and vertical position, velocity, and acceleration traces of the targets. The default settings can be changed by clicking the corresponding cell and choosing a color from the palette or entering the value in RGB or HSB scales;

Acceleration Trace Scaling Factor and **Velocity Trace Scaling Factor**: Set the scaling factor for the acceleration and velocity data so that these two traces can be displayed appropriately in this view;

Message Triangle Width, **Message Triangle Height**, **Button Triangle Width**, and **Button Triangle Height**: Set the width and height of the symbols for message and button events;

Background Color and **Ruler color**: Set the colors for background and ruler of the temporal graph view.

Split Y axis: If true, 0 appear in the middle of the screen, positive values on the top and negative values at the bottom; otherwise, a descending scale is used, with large values on the Y-axis appearing at the bottom of the screen. Note, this setting will only be effective when the data are loaded in a new viewing session.


X Trace Highlighted, **Y Trace Highlighted**, and **Pupil Trace Highlighted**: Whether or not by default the X trace, y trace, and pupil trace are highlighted;

Initial Visible Sample Count: Total number of initially visible samples in the temporal graph view window (the default is 5000).

Use X Position for Fixation Location: If checked, use the X position to indicate the position of a fixation event in the temporal graph view; otherwise, use the Y position.

8.10 Animation Preferences

The parameters related to the animation playback preference settings are covered below:

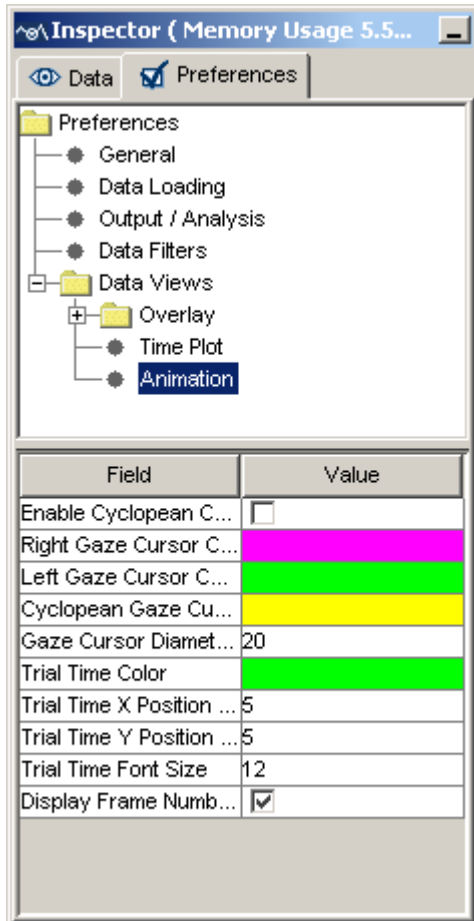

Enable Cyclopean Cursor: For a binocular recording, enabling this option will display a cyclopean gaze cursor (instead of two separate gaze cursors) in the animation view.

Right Gaze Cursor Color, Left Gaze Cursor Color, Cyclopean Gaze Cursor: Set the colors to be used to show the right, left, and cyclopean gaze cursor. The default settings can be changed by clicking the corresponding cell and choosing a color from the palette or entering the value in RGB or HSB scales;

Gaze Cursor Diameter: Sets the diameter of the gaze cursor (in pixels);

Trial Time Color: Sets the color of the trial time code.

Trial Time X Position (%), Trial Time Y Position (%): Sets the x (percent of screen width) and y (percent of screen height) coordinates of the trial time code.

Trial Time Font Size: Sets the font size of the trial time code.

Display Frame Number: If checked, displays the frame number of the video being played.

Display Trial Time: If checked, displays the current playback time relative to the start of the trial.

Display IP Time: If checked, displays the current playback time relative to the start of the current interest period selected.

9 Release History

Version 2.1.1

- General performance improvements and reworked on the playback in the Animation View and video saving.
- Added Dynamic Interest Area Support
 - Modified interest area format
 - Added Mouse Record Mode for creating dynamic interest area instances
 - Added Paste Current and Paste After options
 - Modified interest area creation dialog box
 - Supported position and shape interpolation across instances of Dynamic Interest Area
 - Interest Area data reporting can be either based on the fixation events or samples
- Improved Trial Grouping interface
- Supported adding or removing points for a freehand interest area
- Animation View now displays time relative to start of the trial and/or start of the current interest period.
- Added “Default Video Directory” setting to the General Preferences; supported selection of multiple directories for the “Default Image Directory” and “Default Video Directory”.
- Allowed independent setting for “Hide Spanned Events” and “Trim Spanned Fixation Duration” at the start and end of interest period.
- Added a "Use Substring Match for Messages" option to the Reaction Time and Interest Period definitions.
- Added new variables to the Interest Area Report: IA_DYNAMIC, IA_END_TIME, IA_INSTANCE_COUNT, IA_START_TIME, REPORTING_METHOD, TRIAL_IA_COUNT, TRIAL_TOTAL_VISITED_IA_COUNT.
- Added new variables to the Fixation Report and Saccade Report: CURRENT_/NEXT_/PREVIOUS_FIX_INTEREST_AREA_DATA, CURRENT_/NEXT_/PREVIOUS_FIX_INTEREST_AREA_PIXEL_AREA, CURRENT_/NEXT_/PREVIOUS_FIX_INTEREST_AREA_ID, CURRENT_/NEXT_/PREVIOUS_SACCADE_START_INTEREST_AREA_ID, and CURRENT_/NEXT_/PREVIOUS_SACCADE_END_INTEREST_AREA_ID.
- Added new variables to the Message Report: CURRENT_MSG_INTEREST_AREA_ID.
- Added new variables to the Trial Report: AVERAGE_BLINK_DURATION, and IA_COUNT
- Added new variables to the Sample Report: LEFT_FIX_INDEX, RIGHT_FIX_INDEX, AVERAGE_/LEFT_/RIGHT_INTEREST_AREAS, AVERAGE_/LEFT_/RIGHT_INTEREST_AREA_DATA, AVERAGE_/LEFT_/RIGHT_INTEREST_AREA_DISTANCE,

AVERAGE_/LEFT_/RIGHT_INTEREST_AREA_PIXEL_AREA,
LEFT_SACCADE_INDEX, RIGHT_SACCADE_INDEX

Version 1.11.900:

- Added support for EyeLink 1000 Plus recording files.
- Fixed crashes in loading some EDF files on Mac OSX.
- Added blink correction for pupil size calculation for PUPIL_SIZE_MIN and PUPIL_SIZE_MEAN variables in the trial report; now samples during a blink are not included in the pupil size calculation.
- Added options to exclude some samples before and after a blink event in the pupil size calculation ("Duration to Exclude Before Blinks (msec)" and "Duration to Exclude After Blinks (msec)" in the Data Filters preferences).
- Bug fix for handling blink events in some EDF files.
- Bug fix in handling target position commands ("!V TARGET_POS").
- Supported saving interest area files in UTF-8 encoding.
- Added variables to the Interest Area Report: IA_FIRST_SACCADE_INDEX and IA_LAST_SACCADE_INDEX.
- Bug fix for Interest Area Report variables: IA_FIRST_FIXATION_RUN_INDEX and IA_FIRST_FIXATION_VISITED_IA_COUNT.
- Added option in the "Output/Analysis Preferences" (Identify first/last saccade based on the first last/last fixation) to match the first/last saccade event to the first/last fixation event in the Interest Area Report.

Version 1.11.1:

- Added support for multiple EDF file loading contained in different directories.
- Trial variable value editor now supports copy and paste.
- Added a "Report Hidden" option for the Recording Session and trial groups.
- Added FIXATION_DURATION_MAX, FIXATION_DURATION_MIN, FIXATION_DURATION_MAX_TIME, and FIXATION_DURATION_MIN_TIME to trial report.
- Added IA_LAST_RUN_DWELL_TIME, IA_LAST_RUN_END_TIME, IA_LAST_RUN_FIXATION_%, IA_LAST_RUN_FIXATION_COUNT, IA_LAST_RUN_START_TIME, IA_LAST_FIXATION_RUN, and IA_THIRD_RUN_FIXATION_COUNT to the interest area report.
- Fixed various EDF file loading issues.
- Bug fix for setting the directories in "General Preferences".
- Bug fix for displaying non-ASCII characters in the interest area label, trial variable value editor, and messages.

Version 1.10.123:

- Added support for exporting and importing variable selection in the output reports (trial, fixation, saccade, interest area, sample, and message).

- Added support for exporting and importing interest periods across sessions.
Added support for exporting and importing preference settings.
- Added legend to the grayscale version of fixation maps.

Version 1.10.1:

- This version runs on Windows 2000, XP, Vista, Windows 7, and Mac OSX.
- Better support for non-ASCII characters in the trial variable value editor, interest area label, messages, and report outputs.
- Added "Use UTF-8 Encoding" option to the "Output/Analysis" preference.
- Bug fix for saving the animation view with cyclopean gaze cursor.

Version 1.9.193:

- Added IA_FIRST_FIXATION_X, IA_FIRST_FIXATION_Y, IA_SECOND_FIXATION_X, IA_SECOND_FIXATION_Y, IA_THIRD_FIXATION_X, IA_THIRD_FIXATION_Y, IA_LAST_FIXATION_X, and IA_LAST_FIXATION_Y variables to the interest area report.
- Added CURRENT_MSG_INTEREST_AREAS, CURRENT_MSG_INTEREST_AREA_INDEX, and CURRENT_MSG_INTEREST_AREA_LABEL to the message report.
- Added LAST_BUTTON_RELEASED, LAST_BUTTON_PRESS_TIME, and LAST_BUTTON_RELEASE_TIME to the fixation report. The LAST_BUTTON_TIME variable now returns the time when the last button is pressed or released.
- Added PUPIL_SIZE_MIN_X, PUPIL_SIZE_MIN_Y, PUPIL_SIZE_MIN_TIME, and VISITED_INTEREST_AREA_COUNT to the trial report.
- Bug fix for interest period selection if the end time of the interest period is equal to the start time of a fixation event.

Version 1.9.1:

- Fixation map/heat map is completely redone with the following modifications:
 - Four variants of fixation maps are provided: 1 = based on duration; 2 = based on fixation count; 3 = based on count density; and 4 = based on duration density.
 - A fixation map color legend is provided.
 - Renamed "Minimum Contrast Multiplier" option in the fixation map preference settings to "Low Activity % Cut-off".
 - Supported drawing a title.
 - Transparency setting supported.
 - Option added to customize the Low, Medium, and High activity colors.
 - Option added to set a fixed maximum and minimum value for heatmap scale.

- The compressed EDF files collected on EyeLink CL eye trackers are now supported.
- Added TRIAL_INDEX variable to the fixation, saccade, sample, message, and interest area reports.
- Added IA_FIRST_FIXATION_INDEX, IA_FIRST_FIXATION_RUN_INDEX, and IA_FIRST_FIXATION_VISITED_IA_COUNT variables to the interest area report.