

Research Grade

Isotonic Transducer

User's Manual

Research Grade Isotonic Transducer, 110 VAC/60 Hz 60-3000
Research Grade Isotonic Transducer, 220 VAC/50 Hz 60-3001



HARVARD
A P P A R A T U S

Table of Contents

1

Harvard Apparatus Research Isotonic Transducer User's Manual

SUBJECT	PAGE NO.
Warranty and Repair Information	2
Theory of Operation	3
General Information	4-5
Specifications.....	4
Amplifier/Power Supply	4
Internal Controls and Adjustments	5
Electrical Output	5
Maintenance and Storage	5
Serial Number	5
Operations	6-8
Transducer Setup	6
Smooth Muscle/Isolated Organ Preparation	7
Skeletal Muscle Preparation	8

General Information

2

Serial Numbers

All inquiries concerning our product should refer to the serial number of the unit(s).

Warranty

Harvard Apparatus warranties the instrument(s) for a period of one year from date of purchase. At its option, Harvard Apparatus will repair or replace the unit(s) if it is found to be defective as to workmanship or material.

This warranty does not extend to damage resulting from misuse, neglect or abuse, normal wear and tear, or accident.

This warranty extends only to the original customer purchaser.

IN NO EVENT SHALL HARVARD APPARATUS BE LIABLE FOR INCIDENTAL OR CONSEQUENTIAL DAMAGES. Some states do not allow exclusion or limitation of incidental or consequential damages so the above limitation or exclusion may not apply to you. **THERE ARE NO IMPLIED WARRANTIES OF MERCHANTABILITY, OR FITNESS FOR A PARTICULAR USE, OR OF ANY OTHER NATURE.** Some states do not allow this limitation on an implied warranty, so the above limitation may not apply to you.

If a defect arises within the one-year warranty period, promptly contact **Harvard Apparatus, Inc. 84 October Hill Road, Holliston, Massachusetts 01746-1371** using our U.S. only toll free number 1-800-272-2775 or dial (508) 893-8999. Goods will not be accepted for return unless an RMA (returned materials authorization) number has been issued by our customer service department. The customer is responsible for shipping charges. Please allow a reasonable period of time for completion of repairs, replacement and return. If the unit is replaced, the replacement unit is covered only for the remainder of the original warranty period dating from the purchase of the original device.

This warranty gives you specific rights, and you may also have other rights which vary from state to state.

Repair Facilities and Parts

Harvard Apparatus stocks replacement and repair parts. When ordering, please describe parts as completely as possible, preferably using our part numbers. If practical, enclose a sample or drawing. We offer a complete reconditioning service.



CAUTION: Not for clinical use on human patients.

Theory of Operation

This Isotonic Transducer converts rotary motion into a proportional DC voltage. A moveable capacitor plate attached to the rotating shaft and fixed capacitor plates form a differential capacitor. A patented electronic circuit, (U.S. Patent #4142144) transforms capacitance change linearity into a DC voltage equal to 2.0 volts for 15° of rotation.

The use of the capacitance principle produces the following advantages:

- 1) Extremely low moment of inertia of 0.36 cm which is an order of magnitude better than comparable units. This low inertia insures that complex quick changes are faithfully reproduced.
- 2) Excellent linearity of $\pm 1\%$.
- 3) High voltage output, ± 2.0 volts.
- 4) Excellent stability and freedom from drift.

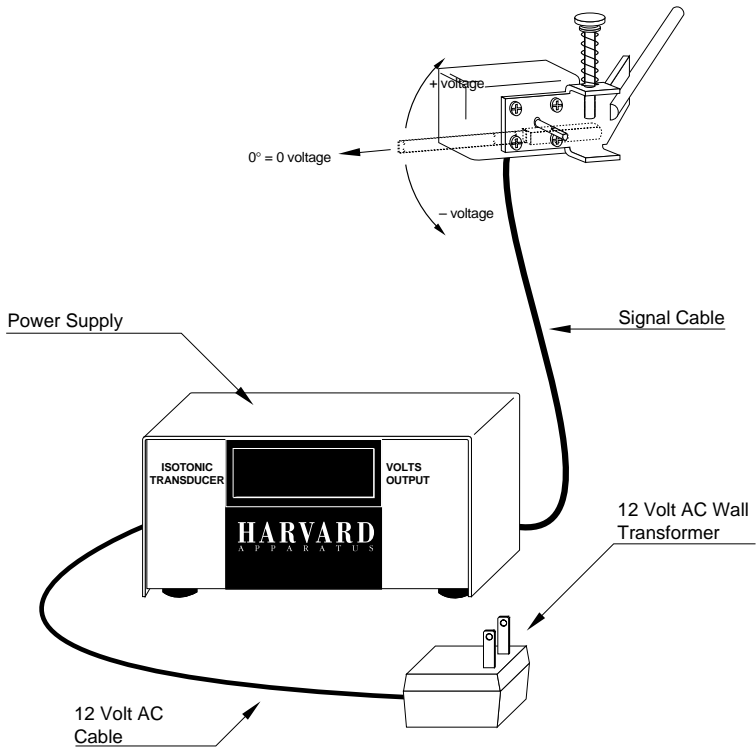


Figure 1. Research Grade Isotonic Transducer

General Information

4

Specifications

Linearity	$\pm 1\%$
Moment of Inertia	0.35 g·cm, without hardware
Breakaway Torque	0.05 g·cm
Accuracy	$\pm 1\%$
Rotation	$\pm 15^\circ$
Output	± 2.0 VDC for $\pm 15^\circ$ rotation
Sensitivity	1 minute of arc
Output Impedance	2 k

Amplifier/Power Supply

- The entire transducer is powered by a 12 volt AC wall transformer.
- There is no ON/OFF switch. The unit can remain powered, as shown on the digital meter, for the entire experiment.

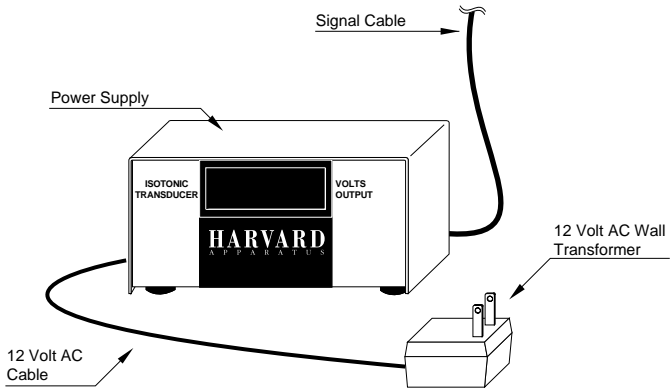


Figure 1. Amplifier/Power Supply

General Information

5

Internal Controls and Adjustments

On the rear panel of the power supply are three small access holes to the trim pots.

The GAIN trim pot is accessed through the center hole and is factory set to deliver 2.0 VDC for 15° of rotation. It can be adjusted to deliver between 1 – 3 volts output for 15° of rotation.

The OFFSET adjustment trim pot is located on the extreme left and has been factory adjusted so that when the flat of the shaft is horizontal, output is zero volts. This can also be user adjusted as required. The right hole is not used.

Electrical Output

- Binding posts are provided on the power supply for output connections. Red is positive (+); black is negative (-) and also ground.
- Output voltage is ± 2.0 VDC for rotation up to $\pm 15^\circ$. Output impedance is 2 k Ω for direct connection to all recorders.

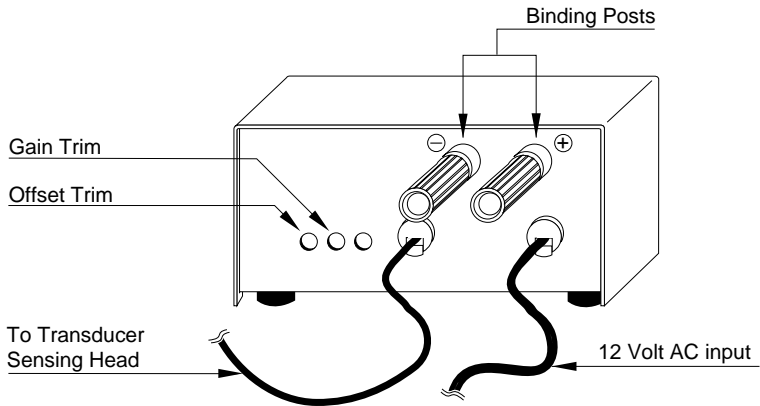


Figure 2. Electrical Output

Maintenance and Storage

- No special precautions are required other than preventing corrosive solutions from entering the space where the output shaft exits the transducer.
- When not in use, store in a clean, dry place.

Serial Number

The serial number is located on the power supply circuit board. Refer to this number in any correspondence.

Transducer Setup

This transducer is supplied with complete hardware to set up the transducer for smooth muscle, isolated organ, or exposed heart applications including bio-assay techniques in an organ bath, or skeletal muscle experiments such as the muscle twitch using the frog gastrocnemius preparation.

The stainless steel mounting rod of the transducer should be securely fastened to a rigid stand for optimal use. It should be noted that maximum sensitivity to muscle movement occurs when the string connected to the muscle is closest to the rotating shaft. In this position, small movements create larger angular motion (see Fig. 1).

When the lever is horizontal and in the same plane as the handle, the electrical output is zero (see Fig. 1). Clockwise rotation produces a positive voltage whereas counter-clockwise rotation produces a negative voltage. The protecting ears of the handle bracket limit rotation to $\pm 18^\circ$. In addition, there are internal limit stops.

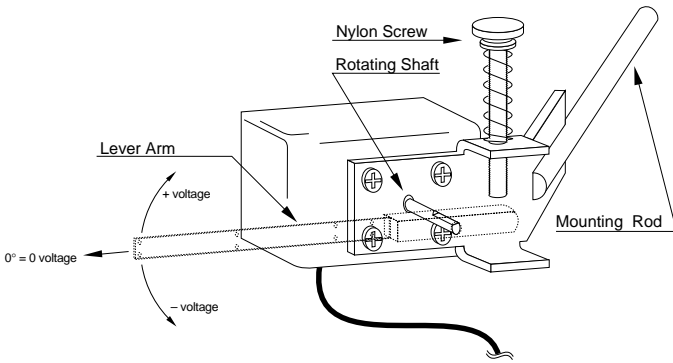


Figure 3. Transducer Setup

Operations

7

Smooth Muscle/Isolated Organ Preparation

For smooth muscle or isolated organs preparations, the following items will be used:

- 1) Simple Wooden or Aluminum Levers
- 2) Lever-Holding Clamp

To use the levers, the lever-holding clamp is attached to the shaft of the transducer using the Allen wrench provided. Make sure that the set screw seats on the flat portion of the shaft. The clamp must be attached to the shaft so that the Allen screw seats on the flattened portion of the shaft. The wooden or aluminum lever is passed through the hole in the clamp and is held in place by the spring of the clamp (see Fig. 2).

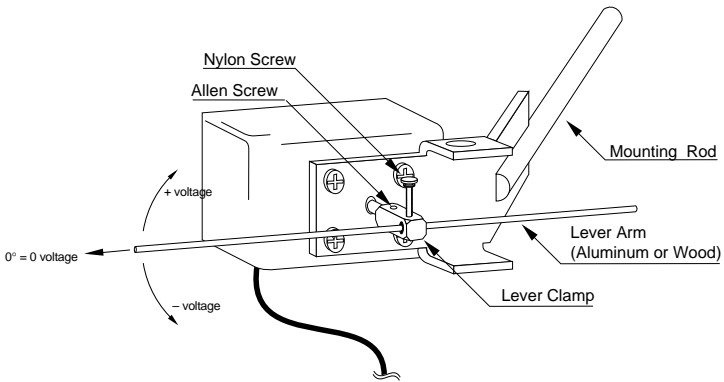


Figure 4. Smooth Muscle/Isolated Organ Preparation

Operations

8

Skeletal Muscle Preparation

For skeletal muscle contraction the following items will be used:

- 1) Muscle Lever
- 2) After Loading Screw and Spring
- 3) Scale Pan
- 4) Weights

Install after loading screw and spring as shown (see Fig. 3). Using an Allen wrench, install muscle lever so that Allen screw bears on shaft. Attach double hook to one of the two positions on the muscle lever and add the scale pan (see Fig. 3).

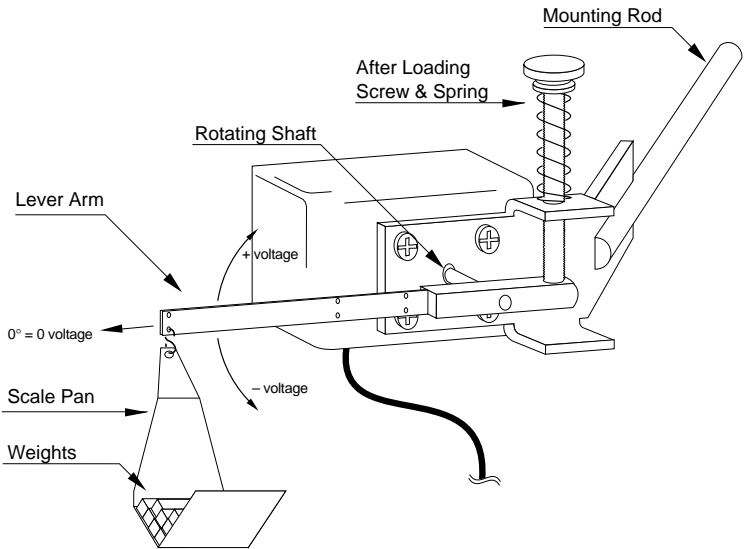


Figure 5. Skeletal Muscle Preparation