



ExpressGene

P5200

Instruction Manual

Release Information

This document is protected by copyright. All Rights Reserved.

Quanta Biotech Ltd

Unit 4 Byfleet Technical Centre

Canada Road

Byfleet

Surrey

KT14 7JX

United Kingdom

Phone +44 (0) 1932 344550 • Fax +44 (0) 1932 353108

Email: info@quantabiotech.com

Web: www.Quantabiotech.com

Manual Version: 1.0

Trademark Information

All product and brand names used in this document are trademarks or registered trademarks of their respective holders.

Copyright © 2010 Quanta Biotech Ltd

TABLE OF CONTENTS

General Information	1
Policy Statement.....	1
Warranty Statement.....	1
1 Safety Warnings and Precautions	2
1.1 What you should know.....	2
1.1.1 Instrument Safety.....	2
1.1.2 Symbols and Conventions	2
1.1.3 Potential Safety Hazard	3
2 Installation.....	6
2.1 Requirements	6
2.1.1 Site	6
2.1.2 Electrical	6
2.1.3 Thermal Block.....	7
2.2 Unpacking the Thermal Cyclor	8
2.2.1 Package Contents.....	8
2.2.2 Initial Inspection	8
2.3 Functional Overview	9
2.3.1 Networking the Thermal Cyclor	10
2.4 AC Power Connection.....	11
2.4.1 Operating Power Requirements	11
2.4.2 Fuses.....	11
2.4.3 Grounding Requirements	11
3 Software Operation.....	13
3.1 Required Computer Skills	13
3.2 Start Screen.....	13
3.3 Administrator Settings.....	15
3.3.1 Cyclor Management.....	15
3.3.2 User Management	17
4 Programming	21
4.1 Creating PCR Programs	21
4.1.1 DND Programming.....	21
4.1.2 Program Wizard.....	27

4.2	General Programming Guidelines	28
4.3	Editing Programs	29
4.3.1	Editing Program Parameters	29
4.3.2	Deleting a Program Step	30
4.3.3	Converting a Gradient Step	30
4.4	Deleting Programs	31
5	PCR Run / Operation.....	32
5.1	User Login	32
5.1.1	Guest Users	32
5.1.2	Registered Users	32
5.2	Changing Users	32
5.3	Selecting and Running a PCR Program	33
5.4	Viewing Script Temperature Profiles Prior to Run	33
5.5	Opening and Closing the Lid	33
5.6	Loading the Cyclor	34
5.7	Setting the Lid Pressure	35
5.8	Program Execution	35
5.8.1	Starting a Run	35
5.8.2	Terminating a Run.....	35
5.9	Reports	37
5.9.1	Viewing Reports.....	37
5.9.2	Deleting Reports	38
5.10	Switching Off the Instrument.....	39
6	USB Memory Stick Operation	40
6.1.1	Importing and Exporting PCR Data Files.....	40
7	Maintenance	41
7.1	Cleaning.....	41
8	Trouble-Shooting	42
8.1	If a Power Failure Occurs.....	42
8.2	Troubleshooting Information.....	43
9	Block Exchange	45

Removing the Block	45
Inserting the Block	45
Appendix.....	46
Instrument Specifications.....	46
Thermal Blocks	46
High Pressure Heated Lid.....	46
Ambient temperature	46
10°C to 30°C.....	46
User Interfaces	47
User & File Management	47
Reports and Validation.....	47
Networking & Barcode Option	48
Thermal Engine Characteristics	48
Power & Dimensions.....	48
Compatible Barcodes.....	49
Compatible Plate Formats.....	50
Order Information	51
Sales & Service	52
Index.	53

General Information

Policy Statement

It is the policy of Quanta Biotech to improve products as new techniques and components become available. Quanta Biotech reserves the right to change specifications at any time.

Warranty Statement

Quanta Biotech guarantees that the thermal cycler you have received has been thoroughly tested and meets its published specification.

This guarantee is valid for 24 months only if the product and functions have been used according to the instruction manual. No liability is accepted for loss or damage arising from the incorrect use of the thermal cycler. Quanta Biotech's liability is limited to the repair or replacement of the unit or refund of the purchase price at Quanta Biotech's option. Quanta Biotech is not liable for any consequential damages.

Quanta Biotech's thermal cyclers are for research use only.

Read the Instruction Manual carefully before using the thermal cycler to ensure that you obtain the best possible results from the machine.

Quanta Biotech thermal cyclers should only be used by suitably qualified and trained people. If the thermal cycler is not used as specified in this Manual, the protection provided by the equipment may be impaired.

1 Safety Warnings and Precautions

1.1 What you should know

Before Operating the Instrument

Ensure that anyone involved with the operation of the instrument is instructed in both general safety practices for laboratories and specific safety practices for the instrument.

Always place the instrument in a location where, if necessary, the main power supply can be disconnected immediately.


1.1.1 Instrument Safety



Please read this section before operating the thermal cycler. Operators of this instrument must be trained in both general laboratory safety practices and the specific safety requirements of the thermal cycler. If the equipment is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.

All functions performed within the context of preparing, performing and completing a run should be done with caution and care, and with general respect both to the instrumentation and to associated chemicals, samples and other devices.

1.1.2 Symbols and Conventions

The following chart is an illustrated glossary of the symbols that are used in this manual.

	CAUTION This symbol indicates a potential risk and alerts you to proceed with caution
---	--

	CAUTION This symbol indicates the presence of high voltage and warns the user to proceed with caution
	CAUTION This symbol indicates risks associated with hot surfaces

1.1.3 Potential Safety Hazard

1.1.3.1 ELECTRICAL

Standard electrical safety precautions should be applied.

- 1 Ensure that the proper voltage is supplied before turning the instrument on for the first time.
- 2 The device must be connected to a grounded socket.
- 3 Do not touch any switches or outlets with wet hands.
- 4 Switch the instrument off before disconnecting the AC power cord.
- 5 Unplug the instrument prior to cleaning up any major liquid spills and prior to servicing any of the electrical or internal components.

Only qualified personnel should perform electrical servicing.

1.1.3.2 EXPLOSIVE SUBSTANCES

Explosive, flammable and reactive substances should never be cycled or incubated in any The thermal cycler.

Do not operate the The thermal cycler in a hazardous or potentially explosive environment.

You must observe the relevant safety regulations when handling pathogenic material, radioactive substances or other substances hazardous to health.

1.1.3.3 PINCHING HAZARD

The lid presents a potential pinching hazard. Do not place hands or fingers into the cycler when the lid is being opened or closed.

1.1.3.4 FLUIDS

Reaction vessels should be filled outside the cycler so that no fluids penetrate the instrument.

Do not submerge the instrument in any liquid.

1.1.3.5 DANGER OF BURNS



Thermal block, inner side of heated lid and reaction vessels quickly attain temperatures of greater than 50°C. **Do not touch – Risk of Burns!** Keep the heated lid closed until temperatures of 30°C or lower are reached.

Do not use any materials (plates, sealings, foils, mats) which are not sufficiently temperature-stable (up to 120 °C).

1.1.3.6 OPERATING ENVIRONMENT

Please ensure that the ventilation slots of the device remain free to vent at all times. A space of at least 10cm should be left around the thermal cycler

The ambient temperature should be between 10°C and 30°C, the humidity between 0% and 95%.

WARNING:

All repair work must be carried out by Quanta Biotech service engineers!

Only original Quanta Biotech replacement parts must be used.

Disconnect the instrument from the mains supply before opening.

2 Installation

2.1 Requirements

The following steps must be followed for optimal and safe operation of the thermal cycler:

2.1.1 Site

- ❖ The thermal cycler should be placed on a rigid, flat, clean surface. Make sure that the instrument stands completely stable.
- ❖ Adequate ventilation is important. Make sure there is sufficient space so that the rear and side air slots are not covered and to allow cooling air to circulate freely around the instrument.
- ❖ There should be no paper under the device as this may block the ventilation path. The unit should always have at least **10cm** distance to the next wall or neighbouring instrument.
- ❖ Quanta Biotech thermal cycler instruments were developed for operation in laboratories in which there is normal ambient temperature and no explosive atmosphere. The ambient temperature should be between 10°C and 30°C the humidity between 0% and 95%.

2.1.2 Electrical

- ❖ The thermal cycler must be connected to a grounded socket.

2.1.3 Thermal Block

- ❖ Reaction vessels should be filled outside the thermal block so that no fluids penetrate the instrument.
- ❖ To prevent damage to the block and the heated lid make sure you use only recommended sample tubes. Unsuitable tubes can be damaged causing the escape of sample material. This may present a serious health hazard, especially when working with infectious materials.
- ❖ Make sure the PCR plates are positioned correctly in the block to prevent damage to the block and the lid.
- ❖ Ensure that the heated lid is closed when heating tubes in the block. Otherwise tubes may burst due to high temperatures thus causing the escape of sample material.
- ❖ The lid presents a potential pinching hazard. Do not place hands or fingers into the cyclor when the lid closes otherwise hands or fingers may be caught between the lid and housing.

2.2 Unpacking the Thermal Cycler

2.2.1 Package Contents

The following parts are included in delivery:

	Q-Cycler II Server
Thermal Cycler	X
Power Cable 110 or 220V	X
Quanta Biotech Thermal Cycler User Manual on CD or memory stick	X

2.2.2 Initial Inspection

All parts of the instrument were thoroughly inspected and tested before the unit was shipped, therefore the instrument should be in good operating order. Carefully inspect the instrument and its accessories for any physical damage sustained in transit.

If the instrument is received in a damaged condition, please file a claim with the transport carrier immediately and contact Quanta Biotech for advice.

If the inspection has passed the Q-Cycler II thermal cycler is ready for use.

2.3 Functional Overview

Quanta Biotech thermal cyclers are automated instruments designed for carrying out the Polymerase Chain Reaction (PCR) and related methods for amplification of DNA templates.



Figure 2-1: the Q-Cycler II (front view)

The Q-Cycler II has a manually operated lid with the ability to adjust pressure for tubes and microplates.

The Q-Cycler II server is operated via a universal software application offering independent control of the server's own thermal block and any combination of up to three, Q-Cycler II, S-96 or S-24 satellite thermal cyclers.



2.3.1 Networking the Thermal Cycler

A network of up to 4 thermal cyclers consisting of one Q-Cycler II server and up to three S-96 and S-24 satellite thermal cyclers can be established or a PC network of up to fifteen S-96 and S-24 satellite thermal cyclers can be established .

USB ports are provided at the side/back of the instruments to connect satellite units to a Q-Cycler II server or a PC via standard USB instrument cables and hub. Optional input devices such as barcode reader, keyboard, mouse or printer can be connected in the same way (see USB connection diagram in Figure 2-4 below).

2.3.1.1 Connecting S-96 and S-24 Satellite Thermal Cyclers to the Q-Cycler II Server or PC.

- 1 Plug the thermal cycler's AC power cord into the mains power connector at the rear of the instrument and connect the other side of the cable with a grounded AC outlet.
- 2 Plug the USB machine cable into the USB port at the side/rear of the Q-Cycler II, S-96 or S-24 satellite thermal cycler.
- 3 Connect the other end of the USB cable directly to a Q-Cycler II server/PC OR to a USB hub, if installing several devices.
- 4 Plug the cable connector of the USB hub into the USB port of the Q-Cycler II server or PC
- 5 Boot up the Q-Cyler Server or PC
- 6 Press the power switch on the instrument(s).
- 7 Launch the application either by powering up the Q-Cycler II server or launching the PCQB application.
- 8 The application will search for Q-Cycler II, S-96 or S-24 satellite thermal cyclers that have been connected and display them in the cycle monitor section of the application. The process can take up to two minutes depending on the number of USB ports in use.

Once found you can configure the thermal cyclers via the **Settings** menu in the navigation bar. For details on software configuration of additional cyclers refer to section Adding a Cycler on page 19.

2.3.1.2 Connecting Optional Devices

Optional input devices such as keyboard, mouse, printer and barcode reader can be connected to the thermal cycler systems via a USB connection.

- 1 Plug the USB cable into the USB port of the input device.
- 2 Connect the other end of the USB cable to:
 - one of the USB ports of the thermal cycler/PC server, **OR**
 - a USB hub, if you wish to connect several input devices to a Q-Cycler II/PC
- 3 When connecting several devices plug the cable connector of the USB hub into the USB port of the thermal cycler server/PC.
- 4 For further configuration of the device consult the manufacturer's instruction manual.

2.4 AC Power Connection

2.4.1 Operating Power Requirements

All Quanta biotech thermal cycler models have self-regulating power. See Instrument Specifications, Power & Dimensions on page 56 for power specifications.

2.4.2 Fuses

WARNING: Disconnect the AC power cord before removing or installing a fuse to avoid the possibility of serious injury from electrical shock!

Fuse compartments are located on the rear of the instrument, above the main power connection. Check the voltage rating label to verify that the instrument is compatible with the AC voltage available at the installation site.

2.4.3 Grounding Requirements

DANGER: please read carefully!

To protect operating personnel, the National Electrical Manufacturers Association (NEMA) recommends that the instrument be correctly grounded. The instrument is equipped with a 3 conductor AC power cord that, when connected to an appropriate AC power outlet, grounds the instrument. To preserve this protection feature, do not operate the instrument from an AC power outlet that has no ground connection.

3 Software Operation

3.1 Required Computer Skills

Users should have basic computing knowledge including using the mouse, selecting items, and drag and drop.

3.2 Start Screen

Shortly after power up of the system, the screen first displays a start-up routine. Then, the Start Screen appears.

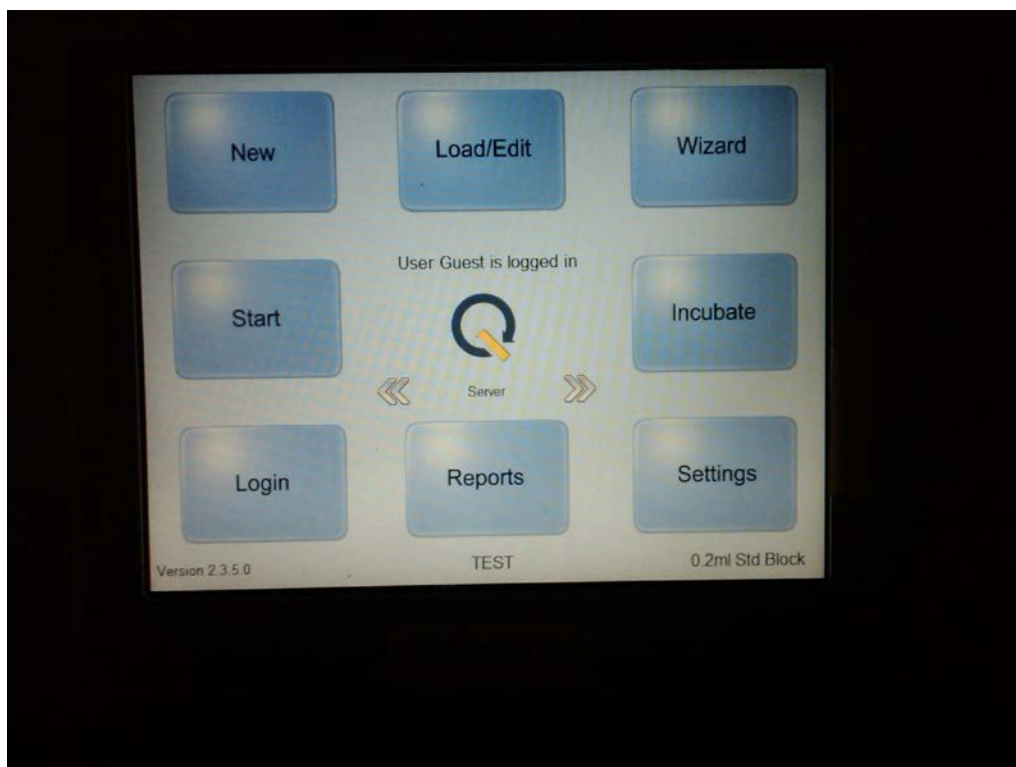


Figure 3-1: Start screen with navigation buttons

Each button describes a task that can be performed upon clicking on it:

New	Creates a New program.
Load/Edit Files	Opens the Explorer to browse stored PCR programs. Use this button to edit or load programs.
Wizard	Opens the Program Wizard dialog window where you can define specific PCR information and a program will be automatically generated
Start	Starts the loaded program. You can view the name of the currently loaded program under the Q logo in the centre of the screen
Incubate	Starts a single temperature incubation for predefined periods or for an indefinite hold.
Login	Enables individual users to login and access there own files and reports
Reports	Opens the report browser where you can search for stored GLP and Lab reports.
Settings	Opens the User Settings dialog window where you can configure user settings such as defining the default start screen and manage cyclers or user data, including power fail and OQ tests.

Q-Cycler II thermal cycler:

The logo in the centre of the screen serves as a return to cycler run button. Clicking on it connects you to the Q-Cycler II run screen for the cycler selected.

The left and right arrows either side of the logo toggle between the server and the different satellites connected.

If the server or satellite name is selected below the Q logo a summary run screen showing all connected blocks appears.

Network Cyclor Monitor					
Cycler Details	Status	Current User	Run Program	Run Information	Time to Completion
Server qCycler	Running	Guest	TEST	Temp: 4.5°C Target: 95.0°C Cycle No: 1	Hours: Min: Sec: 01:46:06
Satellite 2 qCycler Gradient	Running	Guest	TEST	Temp: 4.4°C Target: 95.0°C Cycle No: 1	Hours: Min: Sec: 01:46:06
Satellite 3 qCycler Gradient	Running	Guest	TEST	Temp: 4.9°C Target: 95.0°C Cycle No: 1	Hours: Min: Sec: 01:46:06
Satellite 1 qCycler Gradient	Running	Guest	TEST	Temp: 11.4°C Target: 95.0°C Cycle No: 1	Hours: Min: Sec: 01:46:06

Back

Selecting back takes you to the home screen.

3.3 Administrator Settings

Specific user settings can be defined in the **Settings** menu . Use the **Settings** menu to add new thermal cycler devices or edit existing units, set time and date, or manage user data such as setting up new users, and editing or deleting existing accounts.

3.3.1 Cyclor Management

Q-Cycler II software is designed for use with up to 4 thermal cyclers. Before you can set up additional cyclers in the software, make sure that the new thermal unit is connected to a Q-Cycler II

- 1 To add a cyclor simply connect the new cyclor to the Q-Cyclor II, power up both units and launch the application. The new cyclor will be automatically recognized and installed ready to run. The name and details of the newly connected unit can be changed by an administrator via the settings tab.

3.3.1.1 Editing Cyclor Data

- 1 The name and details of the newly connected unit can be changed by an administrator via the settings button
- 2 Click on **Cyclor Management**
- 3 Click on the thermal cyclor you wish to edit and make the changes
- 4 Confirm your changes by clicking **OK**.

3.3.1.2 Setting Time and Date

- 1 To set the instrument time and date select **Settings**. Then select **Cyclor Settings**.
- 2 Enter the system time and date by pressing the up and down arrow keys.
- 3 Confirm your changes by clicking **OK**.

3.3.2 User Management

Add a new User

User name:

Level:

Password:

Confirm Password:

Virtual Keyboard

1 2 3 4 5 6 7 8 9 0

Q W E R T Y U I O P

A S D F G H J K L

BkSp Z X C V B N M Ok

Space

3.3.2.1 User Levels

With the thermal cycler three different user levels can be distinguished:

- Administrator
- Registered User
- Guest User

Each level has specific user rights. Only administrators are authorized to set up new users and to change and/or assign user rights and passwords.

By default, the thermal cycler starts with Guest user rights.

User Rights	Functions allowed
Guest User	<ul style="list-style-type: none">• Access rights to the Guest folder• Creating, copying, editing and executing programs stored in the Guest folder
Registered User	<ul style="list-style-type: none">• Access rights to the Guest and personal user

	folders <ul style="list-style-type: none"> • Creating, copying, editing and executing programs stored in the Guest and personal folders
Administrator	<ul style="list-style-type: none"> • Access rights to all Guest and user folders • Setup and maintenance of users • Setup and maintenance of cyclers • Archiving/maintenance of all reports

3.3.2.2 First Start-up

Only the administrator is authorized to set up new users and to change or assign user rights and passwords.

As a first-time user you will be logged in as "Guest" user visible in the top status bar. Guest users usually have only restricted user rights and work from the unprotected Guest folder, which can be accessed by all users.

Before you can operate the thermal cycler as a regular user, your administrator has to set you up as a new user. Once you are a registered user, you have a personal, encrypted file folder where you can administer your own PCR program files.

3.3.2.3 Logging in as Administrator

After starting the thermal cycler, the Start screen appears displaying the thermal cycler's navigation elements.

- 1 Click on the **Login** button in the top status bar.
- 2 Select *Administrator* from the **Username** drop-down list and enter the administrator password. The default administrator password is admin
- 1 Click **Login** to confirm your settings.
- 2 You are now logged in as Administrator (visible in the top status bar).

IMPORTANT: It is recommended that the factory-set login password be changed to prevent unauthorized access to the instrument

3.3.2.4 Adding or Modifying User Accounts

NOTE: Only the administrator can set up new users or modify existing user accounts.

Adding New User Accounts

- 1 To add a new user, click **Settings**. The **User Settings** dialog opens.
- 3 In the **User Management** section click **OK**.
The **User Management** dialog opens listing the available user names and corresponding user levels.
- 4 Click **New User** to open the **Add a new User** dialog.
- 5 Assign a user name and password..
- 6 Select the required user level (Guest, User, Administrator) from the drop-down list.
- 7 Now you can define the user's script reading and writing privileges by checking the appropriate boxes:

Edit/Rename Script	The user has the right to edit or rename a program file
Delete Script	The user has the right to delete program files
Create Script	The user has the right to create program files
- 8 Once you have entered all data, click on **Create User** to save the settings. After successful setup, a confirmation message will be displayed and the new user name will be added in the **User Management** list.

Modifying User Accounts

- 1 To modify an existing user account click **Settings**. The **User Settings** dialog opens.
- 9 In the **User Management** section click **OK**.
The **User Management** dialog opens listing the available user names and corresponding user levels.
- 10 Click **Edit User** to open the **Edit a User** dialog.
- 2 Here you can modify user data such as name, user level, password, user privileges and script directory protection.

- 3 Click **OK** to save the changes.
A confirmation message will be displayed after successful update of the user data.

Deleting User Accounts

NOTE: User accounts can only be deleted by the administrator.

- 1 To delete an existing user account click **Settings**. The **User Settings** dialog opens.
- 2 In the **User Management** section click **OK**.
The **User Management** dialog opens listing the available user names and corresponding user levels.
- 3 Select the user name you wish to remove.
- 4 Click on **Remove User** to remove the user name from the user list. It removes all user permissions that the user has and does not allow that user to access any the thermal cycler device. It also deletes all scripts and GLP reports associated with that user.

NOTE: The Remove User command also deletes all scripts and GLP reports associated with that user. The Administrator user account cannot be deleted.

4 Programming

4.1 Creating PCR Programs


4.1.1 DND Programming

Drop-and-Drag programming (DND programming) is an easy and intuitive method for conveniently creating PCR programs step-by-step.

The basic idea of DND programming is that you simply choose the required command from a command list, drag it with the mouse into the program window on the left and drop it into the desired program position. This opens a dialog window where you can define further command-specific settings such as time, temperature, lid pressure, etc. In this way you can create your PCR program command by command according to the planned programming sequence.

The command order can be changed easily any time just by dragging the selected command on the desired position within the program window.

Once the program is started the cycler will process the commands from top to bottom according to the defined programming sequence.

- 1 To generate a new PCR program using DND, just click  on the Start screen. This will open the **Compose a New PCR Program** dialog with its drag-and-drop functions:



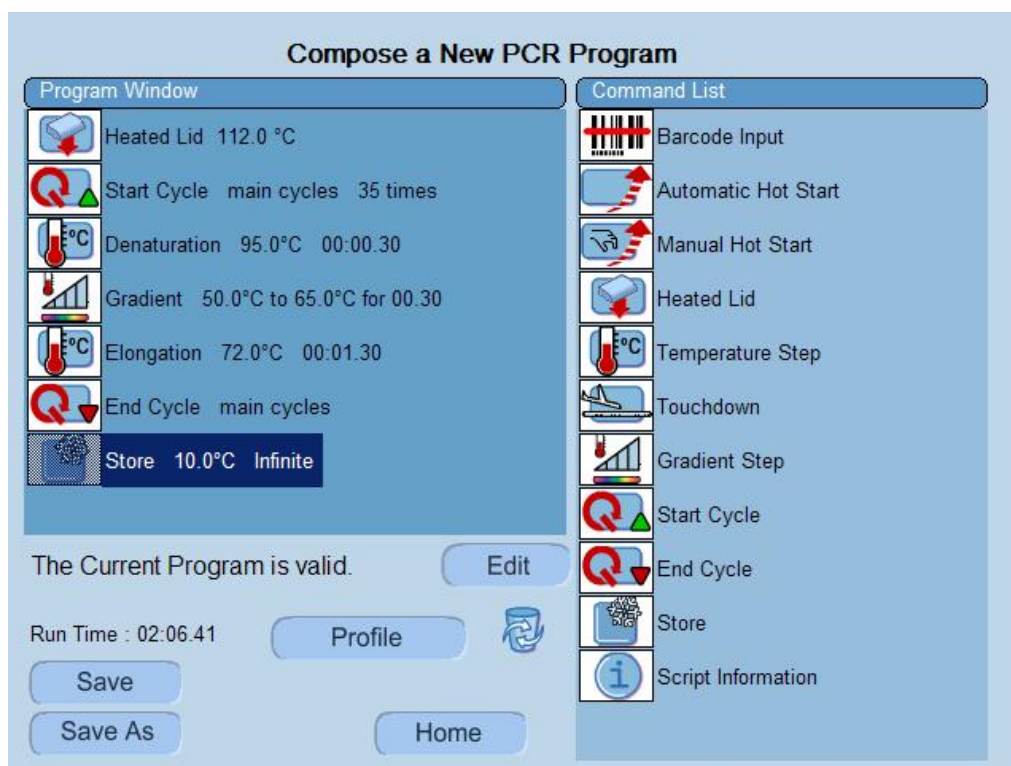


Figure 4-1: Compose a New PCR Program dialog

The dialog window consists of the *Command List*, which contains all available program steps, and the *Program Window* where you can create or edit PCR programs.

- 11 Select a command from the *Command List*, then drag and drop it into the *Program Window*. This will open the dialog window of the selected command.
- 12 Enter your settings in the dialog fields of the selected command.
- 13 Confirm your settings by clicking **OK**. This will return you to the **Compose a New PCR Program** window.
- 14 To add another command to your program repeat steps 2–4 for each required command.
- 15 Once you have finished creating your program, save it by clicking on **Save As** and you will be prompted for a file name.
- 16 Name your file whatever you like within Windows naming conventions.
- 17 Click on **Profile**, if you wish to view a graphical representation of the defined temperature profile.

4.1.1.1 Program Steps

The following program steps are available in the Command List for creating customized PCR programs in the DND environment:



Barcode Input

Use this command to define barcode settings. When composing the program, it is recommended to place the barcode command at the beginning. During program run you will be asked to read the plate's barcode. You may do this with a handheld USB barcode scanner or by manually entering the barcode. If you use the barcode scanner, the barcode cannot be edited after reading.

The program will continue after you have placed the plate into the cycler and pressed *OK*. The barcode will be recorded in the GLP and LabBook report and displayed in the *Info* section of the Cycler screen.

Scan Using Barcode Entry: Barcode reading using the barcode scanner

Scan Using Keyboard Entry: Manual entry of the barcode via keyboard

Path for worklist: Specifies the path for the worklist.



Automatic Hot Start

Automatic Hot Start takes advantage of novel Polymerase technology, which enables the performance of an extended hot start without degrading the enzyme at high temperature. This is realized e.g. using a bead or antibody protection. The protection is released after the extended denaturation is elapsed and the enzyme starts working.

When using the *Automatic Hot Start*, you just have to select the denaturation temperature and time. Running the program, the next command will be executed immediately after the denaturation time is elapsed.

Denaturation Temperature: Specified denaturation temperature

Hold for: Specified time of the program step

Use with automatic Hot Start Polymerase: Examples for denaturation times

Confirmation Change Polymerase 900 sec denaturation

Please also refer to the polymerase manufacturer's recommendations



Manual Hot Start

When selecting the *Manual Hot Start* command, you just have to supply the denaturation temperature and time. The instrument will notify you after the denaturation time is elapsed. You will then be asked to open the cycler and add the enzyme. The instrument timer is now on hold and will continue only after you have pressed *OK* in the displayed *Hot Start* information window.

Denaturation Temperature: Defined denaturation temperature

Hold for: Specified time of the program step

Lid

Use this command to set the required lid temperature.

Lid Temperature: To determine the temperature of the heated lid. The temperature may range from 100°C to 112°C. The minimum increment is 1°C.

Lid is ON: Select this option to activate lid heating.



Temperature Step

Use this command to increment the processing temperature (in °C) by a preset value and to define the step duration.

Processing Temperature: Specified processing temperature in °C.

Hold for: Specified time of the temperature step.

Temperature Ramp: Rate of change of the sample temperature. Rates can be from 0.1°C to 3°C per second. When setting the ramp to zero, the instrument ramps with maximum capacity.

Temperature Increment: Temperature increment and decrement per cycle. Steps can be from 0.1 °C to 10 °C.

NOTE: For decrement use negative figures (e.g. -2)!

Time Increment: Time increment per cycle in seconds.

Step Type: Indicates the selected PCR step type for information purposes. The supplied step type information will then be displayed in the online cycler monitor.



Touchdown

Touchdown is an advanced PCR technique used to reduce unspecific binding. With the thermal cycler, Touchdown PCR is provided as a one-step command for easy implementation in the programming sequence.

Denaturation Temperature: Defined denaturation temperature.

Maximum Annealing Temperature: Sets the maximum annealing temperature.

Minimum Annealing Temperature: Sets the minimum annealing temperature.

Hint: The temperature difference is commonly between 5°C and 10°C starting with a temperature that is 2°C above the higher Primer melting temperature. Example: if T_m Primer 1 is 60°C and T_m Primer 2 is 54°C, the resulting Touchdown will be from 62°C to 52°C.

Elongation Temperature: Sets the elongation temperature.

Cycles: Number of required cycles for the Touchdown section.

Duration: Duration of the temperature step in seconds.



Gradient Step

Use this command to define settings for the gradient function. Just enter the required minimum and maximum temperature gradient and time.

Minimum Temperature Gradient: To set the minimum temperature of the gradient step. Min. temperature must not be below 30°C.

Maximum Temperature Gradient: To set the maximum temperature of the gradient step. Max. temperature must not exceed 80°C.

NOTE: The maximum / minimum gradient span is 20°C / 4°C.

Hold for: Duration of the Gradient step.

Convert Gradient Step to a standard temperature step: The Gradient step can be converted easily to a temperature step after the annealing temperature optimization experiment. To do so, just check the checkbox and select the temperature column with the best result of your optimization experiment. The converted gradient step will then automatically use the temperature of this best-performing column.

Gradient steps are not available for the S-24 thermal cycler

NOTE: Resulting temperature steps may be reconverted to gradient steps any time by disabling the *Conversion* option



Start Cycle

Starts a PCR cycle.

Cycle Name: Enter the name of the cycle.

Number of Loops: Enter the required number of loops.



End Cycle

Insert this command to terminate a PCR cycle.



Store

Cools the thermal cycler down to a temperature between 4°-12°C for a specified period of time.

Store Temperature: Storage temperature in °C.

Hold for: Duration of the Store step.

Infinite: Holds the specified temperature for an unlimited period of time.

NOTE: When executing the *Store* command, the lid pressure is set to zero Newton



Script Information

Use this command to record your comments about the PCR program.

4.1.2 Program Wizard

The Program Wizard™ is a unique feature that makes it easy for first-time PCR users to write their own protocol. The Program Wizard takes care of the calculations required for generating a PCR program. All you need is to supply the biological data of the Primer and template DNA such as Primer sequence or melting temperature, template length and origin of DNA.

The Program Wizard is also an ideal tool for first shot PCRs as it uses proven calculations, e.g. for the Annealing Temperature.

- 1 To generate a new PCR program using biological data, click on **Program Wizard** in the start screen.
- 2 The **Primer Information** dialog opens from where you will be guided step by step through the programming procedure.
- 3 After you have supplied all required data click on *Generate PCR Program*. The **Compose a New PCR Program** dialog window opens where you can view and edit the previously defined program steps.
- 4 To edit a program step, double-click the respective command.
- 5 To view a graphical representation of the defined temperature profile, click **Profile**.

4.2 General Programming Guidelines

To ensure a correct PCR run, the following guidelines should be followed when composing a new program:

- A program must contain at least one *Temperature Step*.
- Cycles cannot be nested. A started cycle must have an *End Cycle*

command before starting the next *Start Cycle* command.

- *Barcode Input* command cannot be programmed inside a cycle.
- Only a single *Script Information* command is valid within a program.

The following commands or command combinations are considered "stages":

- ▶ Automatic Hot Start
- ▶ Manual Hot Start
- ▶ A string of up to ten single Temperature steps
- ▶ Single Gradient steps
- ▶ A complete cycle including Temperature or Gradient steps within a start cycle / end cycle pair.
- ▶ A Touchdown command
- ▶ A Store command

When using these stages in a script, the following rules apply:

- The maximum number of stages within a script must not exceed 10.
- The maximum number of Temperature and Gradient commands inside a cycle must not exceed 10.
- The maximum number of cycles/loops within a stage is 99.

4.3 Editing Programs


4.3.1 Editing Program Parameters

You can edit the parameters of existing programs and store the program with the same name or rename the edited program.

To edit an existing PCR program click on the load/edit button on the main screen:

- 1 Depending on your user level (guest or registered user), the specific user directory or guest directory is displayed listing all existing PCR program files.
- 2 Select the program you wish to edit.
- 3 Click **Open**.
Opening the file automatically loads the PCR program and the program steps are displayed in the Program Window of the **Compose a new PCR Program** dialog.
- 4 Double-click the command you wish to edit. This will open the dialog window of the respective command.
- 5 Edit the parameters as needed.
- 6 Click **Save** to save your changes and to return to the **Compose a new PCR Program** window.
- 7 Repeat steps 4–6 for each command you wish to edit.
- 8 After you have modified all data, click **Save** in the **Compose a new PCR Program** dialog, if you wish to store your changes in the existing PCR program. OR
- 9 Click **Save As**, if you wish to rename the program. You will then be prompted for a file name.

4.3.2 Deleting a Program Step

- 1 To remove a program step from the program just drag it from the Program Window and drop it into the recycle bin  icon located below the Program Window.
- 2 Save the program as described above.

4.3.3 Converting a Gradient Step

A gradient step can be converted to a standard temperature step with a few mouse clicks.

- 1 After performing the gradient experiment, have a look at the gel image and select the best-performing column.
- 2 Double-click the Gradient Step command to open its dialog window.
- 3 Select the **Convert Gradient Step to a standard temperature step** checkbox.
- 4 Select the appropriate column from the drop-down list in the Conversion section.
the thermal cycler software will now convert the gradient step into a temperature step and automatically select the right temperature in the chosen column.

NOTE: If it is required to repeat the gradient experiment, the converted temperature step can be set back to a gradient command just by unchecking the Conversion option

4.4 Deleting Programs

- 1 To delete an existing PCR program Click Load Edit in the start screen.
- 2 Depending on your user level (guest or registered user), the specific user directory or guest directory is displayed listing all existing PCR program files.
- 3 Select the program you wish to remove.
- 4 Click **Delete**. This will remove the program from the script folder.

5 PCR Run / Operation

5.1 User Login

5.1.1 Guest Users

When powering up the instrument the default user level is Guest. If you wish to continue to operate the cycler with Guest user rights, no special login procedure is required. In this case you can access the unrestricted Guest folder and operate the cycler as described in the relevant chapters.

5.1.2 Registered Users

Once the administrator has set you up as a registered user (see 3.3.2.4), you can operate the thermal cycler using your personal data and file folders.

To log into the system as a registered user proceed as follows:

- 1 Click on the **Login** button in the main screen– the **Login** dialog opens.
- 2 Select your user name from the Username drop-down list and enter your password.
- 3 Click on the **Login** button. This will log you into the system giving you reading and writing rights as specified for your user level.
- 4 After successful login your user name is displayed in the top status bar.

NOTE: In case a user forgets his or her password, the logged-on administrator can assign him/her a new one.

5.2 Changing Users

In order to log into the system when another user is already logged in, proceed as described in section *Registered Users* above.

5.3 Selecting and Running a PCR Program

- To select an existing PCR program choose **Start** from the main screen
- 1 Your personal file folder opens or, if you are logged in as a guest, the guest folder opens displaying the available PCR programs.
 - 2 Select the program you wish to run.
 - 3 Click **Run selected** on the right-hand side of the window to start the selected program.

NOTE: The *Start PCR* command button in the navigation bar on the left-hand side of the screen will start the currently loaded program, e.g. the previously used program, but not the selected program!

- 4 If you wish to view the PCR program steps, click **Open**.
Opening the file automatically loads the PCR program. The name of the loaded program is displayed in the top status bar.
- 5 After a program is loaded you may view its script parameters prior to run by clicking on **Profile**.

5.4 Viewing Script Temperature Profiles Prior to Run

Before starting a run, you can view the temperature profile of a new or stored program.

- 1 If not yet displayed, open the desired program as described in section 5.3, *Selecting and Running a PCR Program*.
- 2 Click on **Profile** in the **Compose a New PCR Program** dialog window.
An information screen is displayed detailing all the parameters of the selected program.

5.5 Opening and Closing the Lid

For Q-Cycler II thermal cyclers the lid can be opened and closed manually at any stage.

5.6 Loading the Cyclor

For Q-Cyclor II thermal cyclers open the lid manually.

WARNING:



Danger of Burns! The thermal block, test tubes and plates may reach temperatures as high as 100°C! Before loading or unloading the cyclor, keep hands away until temperature reaches 30°C or less

Pinching Hazard! The lid presents a potential pinching hazard

D o not place hands or fingers in to the cyclor when the lid closes

Loading Plates

Depending upon the thermal block installed, the thermal cyclor can be loaded with one 96-well PCR plate or one 384-well PCR plate.

Put your plate on the block and ensure that it is correctly positioned in the provided space.

Loading Tubes

Q-Cyclor II thermal cyclers can be loaded with up to 96 PCR tubes (0.2 ml) or 48 PCR tubes (0.5 ml) when using the 96-well combi block.

Distribute tubes equally over the block.

NOTE: When using a Gradient step inside your program, use the central rows in the block (rows D + E).

5.7 Setting the Lid Pressure

Simply close the lid and turn the knob. This sets the pressure to low pressure for tubes. If high pressure for plates is required simply turn the knob slightly further

5.8 Program Execution

5.8.1 Starting a Run

To start a run click start from the main screen

Your personal file folder opens or, if you are logged in as a guest, the guest folder opens displaying all available PCR programs.

- 1 Select the program you wish to run.
- 2 Click **Run selected** on the right-hand side of the window to start the selected program.

NOTE: The *Start PCR* command button in the navigation bar on the left-hand side of the screen will start the currently loaded program, e.g. the previously used program, but not the selected program!

After a run is started, a graphical representation of the temperature parameters is displayed. In addition, detailed program and instrument information can be viewed using the **Program** or **Info** buttons located in the right-upper corner of the information screen.

Click **Program** to display the program steps of the running protocol. The executed step is highlighted.

Click **Info** to view detailed instrument information such as temperature sensor values, lid status, remaining time and step time or GLP file name.

5.8.2 Terminating a Run

After a run is started, the **Start PCR** button in the start screen change to **Stop PCR**.

- 1 To stop a program manually before it completes, ensure the appropriate thermal cycler is selected and press **Stop PCR**.

To open the cycler turn the knob located on the top of the instrument
The manual abortion of the program will be recorded in the GLP and lab book reports.

5.9 Reports

Browse Generated Reports [Archive Reports](#)

☒ Administrator
☒ Guest

Report Name	Creation Date
Incubation.glp	01/October/2007 at 1
Incubation.lab	01/October/2007 at 1
pcr optimisation.glp	01/October/2007 at 0
pcr optimisation.lab	01/October/2007 at 0
TST1.glp	28/September/2007 a
TST1.lab	28/September/2007 a
Incubation.glp	27/September/2007 a
Incubation.lab	27/September/2007 a
Incubation.glp	27/September/2007 a
Incubation.lab	27/September/2007 a
TST1.glp	27/September/2007 a
TST1.lab	27/September/2007 a
Incubation.glp	27/September/2007 a
Incubation.lab	27/September/2007 a

☒ Group by Date
☐ Group by Program
☐ Group by Cycler

☒ Show GLP ☒ Show LabBooks

[View](#)
[Home](#)

GLP Report
c:\qCycler\Users\Guest\GLP\TST128092007133700701.glp
Barcode: None

Date 28/09/2007 Run Time 00 days 01:22:39 Program TST1
User Guest Cycler qCycler
Last Calibration 27/09/2007 Next Calibration 26/11/2007 Serial 701

Run was terminated Manually

Step	Stage	Step	Cycle	Temperature	Duration	Time
TD Denaturation 95.0°C ...	1	1	2	94.9	00:30	0:06.0
TD Annealing 57.9°C 00:30	1	2	2	58.1	00:30	0:06.4
TD Elongation 68.0°C 0...	1	3	2	68.0	04:00	0:10.4
TD Denaturation 95.0°C ...	1	1	3	94.9	00:30	0:11.3
TD Annealing 57.4°C 00:30	1	2	3	57.4	00:30	0:12.2
TD Elongation 68.0°C 0...	1	3	3	68.0	04:00	0:16.3
TD Denaturation 95.0°C ...	1	1	4	95.0	00:30	0:17.3

[Back](#)

5.9.1 Viewing Reports

Depending on the assigned user level, you are allowed to view and print LabBook reports stored in the following file folders:

User rights	File folder	Action
Guest users	Guest folder	Viewing and printing Lab Book reports The access to GLP reports is denied .
Registered users	Guest folder & Personal file folder	Viewing and printing Lab Book reports The access to GLP reports is denied
Administrator	All guest and user folders	Viewing Lab Book reports Access to GLP reports permitted

To view reports proceed as follows:

- 1 Click **Reports** in the navigation bar. This will open the **Browse Generated Reports** window listing the generated reports.

- 2 Group the displayed reports by selecting one of the sorting criteria below:

Group by Date: Sorts records in ascending order by date.

Group by Cyclor: Sorts records by cyclor name.

Group by Program: Sorts records by program name.

- 3 For viewing GLP and/or LabBook reports select one of the following options:

Show GLP: Displays the generated GLP records (only for users with administrator rights)

Show LabBooks: Displays the generated LabBooks

5.9.2 Deleting Reports

To delete reports proceed as follows:

- 1 Open the **Reports** dialog window as described above.
- 2 Select the report you wish to remove.
- 3 Click **Delete** to remove the report from the reports list.

NOTE: Only users with administrator rights are authorized to delete reports.

5.10 Switching Off the Instrument

The unit can be switched off at any time during operation using the power button at the front of the instrument. There is no need to exit the thermal cycler software prior to switching off the instrument.

6 USB Memory Stick Operation

6.1.1 Importing and Exporting PCR Data Files

You can import Primer sequences and PCR programs created on a different thermal cycler or on a PC, or export PCR programs and reports that have been created using a USB memory stick.

Exporting Program Files to a USB Archive

- 1 Make sure the USB memory stick is inserted in the USB slot on the front of the instrument.
- 2 Click **View & Edit Files** to display your personal file folder. Select the desired PCR program as described in section 5.4 Selecting a PCR program.
- 3 Click **Save As** to operate the **Save Program** dialog window.
- 4 Enter the name of the program.
- 5 Activate the **Save to USB** checkbox. The file is then saved to the USB archive.

NOTE: the lock button on the memory stick may prevent data access. Please ensure the lock button is in the unlocked position.

Importing Files from a USB Archive to the Thermal Cycler

- 1 Make sure the USB memory stick is inserted in the USB slot on the front of the instrument.
- 2 Click **View & Edit Files** to display your personal file folder. Click **View USB** to access the USB archive.
- 3 Select the desired file from the archive.
- 4 Click **Save As** to select a file name and path.
- 5 Confirm with **OK** to save the file.

7 Maintenance

7.1 Cleaning

Quanta Biotech thermal cyclers are designed to require a minimum amount of maintenance by the user. They can be cleaned using water or a mild laboratory-cleaning agent.

The instrument should not come into contact with organic solvents or aggressive solutions. Ensure that no liquid enters the thermal cycler. For safety reasons, the device must be switched off and disconnected from the power supply prior to cleaning.

The electrical safety fuses are located above the main power switch and the main power plug at the rear of the device. Before the fuses are replaced, the device must be switched off and disconnected from the mains supply. Only fuses with the correct voltage values may be used.

Only qualified service personnel should perform servicing of the thermal cycler.

No warranty obligation is assumed by Quanta Biotech in the event of damage caused by unauthorized servicing.

8 Trouble-Shooting

8.1 If a Power Failure Occurs

All Quanta Biotech thermal cyclers are equipped with an autopower fail restart function that allows for power failures and safe continuation of a PCR run after resumption of power if set to do so in the cycler settings.

The autopower fail restart function restarts the script from the beginning of the temperature step in which it failed.

The following table lists the three power failure scenarios and the actions the instrument takes if the power is interrupted during operation.

Observation	Explanation / Remedy / Actions
The computer fails but the thermal cycler unit still has power	In this case the thermal engine will continue to execute the script without any interruption.
The thermal cycler fails but the computer still has power	The execution of the script will be affected and will halt whilst the power is off. Once the power resumes, restart the unit and the script will resume from the beginning of the step it was executing when the power failed.
Both the computer and the thermal cycler unit lose power	<p>The execution of the script will be affected and will halt whilst the power is off. Once the power resumes the script will resume from the beginning of the step it was executing when the power failed.</p> <p>Corrective action:</p> <p>Restart the application and the connection will be restored.</p>

8.2 Troubleshooting Information

Refer to the following table for a description of potential problems and recommended actions that you should take.

Observation	Explanation / Remedy / Actions
No actual temperature displayed	Block not correctly inserted in the machine. Corrective action: Check that the block is inserted correctly and reboot the machine.
One thermal engine thermistor does not agree with the others	Run Thermal Engineer OQ test ⁽¹⁾ to assess whether performance is within specified limits.
Poor experimental yield	1. Run Thermal Engine OQ test test ⁽¹⁾ to check that thermal engine operation is within specification. 2. If transferring protocol from a different make of thermal cycler, use TAS ⁽²⁾ system to determine new protocol parameters.
Program does not start	Block not correctly inserted in the machine. Corrective action: Check that the block is inserted correctly and reboot the machine.
Tubes are crushed	Applied pressure too high. Corrective action: Reduce pressure applied using lid knob
Condensation in tubes during a run	Condensation will appear naturally at the end of a run and has no deleterious effects. Simply spin or flick the liquid back to the bottom of the tube and continue with the post PCR steps. However, condensation during a run will result in sub-optimal cycling. If this is experienced try increasing the lid temperature.
Sealed microplate losing sample volume	Poor quality microplate/sealing system or inadequate pressure. Corrective action: Increase the pressure by turning the knob in the lid
USB devices including memory stock and barcode scanner not recognised	Reboot machine.
Satellites are not recognized by the application	Check all satellites are powered and connected. reboot Server.

Observation	Explanation / Remedy / Actions
Machine does not power up	Check power is on at supply. Check fuses at rear of instrument above power inlet.
Blank Display lit but showing no information	Reboot machine. If the problem persists there is an internal connection problem: call Technical Support.
Abnormal block heating/cooling performance	Run Thermal Engine OQ test ⁽¹⁾ to check performance.

(1) Please refer to the thermal cycler IQ/OQ Protocol document for instructions on performing instrument qualification tests.

(2) The Temperature Acquisition System (TAS) can be purchased separately from Quanta Biotech Ltd.

9 Block Exchange

Block exchange can be performed within seconds. No tools are required for this purpose.

CAUTION:



Turn power off and interrupt power line before changing the block



Removing the Block

Remove the block by simply holding it with two hands on the left and right side and lift it up.

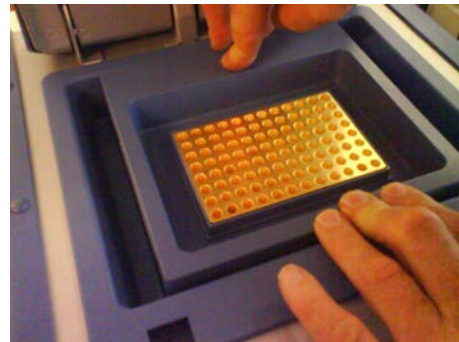
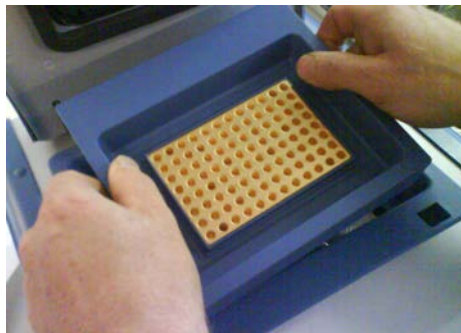


Figure 9-1: Exchanging the block

Inserting the Block

Hold the block with two hands on the left and right side, with the block connector at the back of the block, and place it into position in the thermal cycler.

Appendix

Instrument Specifications

Thermal Blocks

Block Materials	Modular anodized aluminium or gold coated silver blocks with 4 sensors
Traceability	NIST traceable temperature calibration
Blocks Available	<ul style="list-style-type: none">- 96 well gradient block: For 96 well PCR plates or 96x 0.2 ml tubes- 96 well gradient combi block: For 96 well PCR plates, 96 x 0.2 ml tubes- 4 slide/microarray- 384 well high throughput block: for 384 well plates

High Pressure Heated Lid

Lid Temperature Range	To 112°C, user selectable in 1°C increments
Lid Pressure Range	Lid pressure manual - Compatible with all seals
Ambient temperature	10°C to 30°C
Relative Humidity	0% to 95%

User Interfaces

Screen Type	6.4 inch TFT colour touch-screen
System Input	Touchscreen, external keyboard & mouse (USB), barcode scanner (USB)
Temperature Display	Real-time graphical display of actual & set block temperatures Real-time digital display of all sensor values
Communication Interfaces	1x USB front (personal memory stick)

User & File Management

User Level Management	3 levels: Administrator, User with administrator selectable restricted rights, Guest
User File Protection	Yes, for programs and reports
File Organization	Windows® Explorer, user-defined folders & subfolders
Program Storage	Minimum 1000 (internal memory or USB stick). Unlimited PC program storage

Reports and Validation

Report Function	Encrypted GLP report, LabBook report
Validation	Internal auto validation prior to each program start Thermal engine validation function

Networking & Barcode Option

Networking	Cycler network up to 4 units Server - Satellite system network up to 4 units
Barcode Option	Barcode documentation via handheld barcode scanner

Thermal Engine Characteristics

Temperature Control	4°C - 99.9°C with simulated volume dependent control algorithm
Sample Volume Range	5-100µl
Sample Accuracy	+/- 0.4°C (20-99°C), +/- 1°C (4-20°C)
Sample Homogeneity	+/- 0.4°C
Heating	Up to 5°C per second
Cooling	Up to 3°C per second
Sample Overshoot	< 1°C
Gradient Temperature Range	30°C - 80°C (96 well only)
Maximum/minimum Gradient Span	30°C / 4°C

Power & Dimensions

Electronic Power Supply	100V to 240V (frequency 48 to 62Hz)
Dimensions	(WxDxH) 34cm x 42cm x 26cm
Weight	12 kg approx
Operating conditions	10°C - 30°C, 0 - 95% relative humidity
Regulatory	CE, UL, CSA compliant
Warranty	2 years on all systems

Compatible Barcodes

The following barcodes are compatible with THE Q-Cycler II's documentation system:

- 2 of 5 interleaved
- Codabar
- Code 128
- Code 39
- Code 39 full ASCII
- Code 93 full ASCII
- EAN-13, EAN -13 with 2 supplement, EAN -13 with 5 supplement
- EAN-8, EAN -8 with 2 supplement, EAN -8 with 5 supplement
- EAN-128
- UPA, UPA -13 with 2 supplement, UPA -13 with 5 supplement
- UPE, UPE -13 with 2 supplement, UPE -13 with 5 supplement
- MSI
- C.I.P

Compatible Plate Formats

The following plate formats are compatible with the thermal cycler:

Manufacturer	Art. No. & Product Name	Type
ABgene		
	AB 800	96 well, skirted
	AB 900	96 well, semi-skirted
	AB 1111	384 well
	TF-0384	384 well
Bilatec		
	THERMOSPRINT® 96 PP, transparent, 311296; only w/o workstation	96 well, non-skirted
	THERMOSPRINT® 384 PP, transparent, 310384	384 well
NUNC		
	Art. No.: 240600, 96 well, non-skirted.	96
	Art. No.: 230013, 96 well, semi-skirted	96
	Art. No.: 230012, 96 well, skirted	96
	Art. No.: 230584, 384 well	384
Eppendorf		
	Art. No.: 0030 128.648, twin tec PCR plate, 96 skirted	
	Art. No.: 0030 128.575, twin tec PCR plate, 96 semi-skirted	
	Art. No.: 0030 128.508, twin tec PCR plate, 384	
Corning		
	Thermowell® 96 Well Polypropylene PCR Microplate, Natural (#6551)	
	Thermowell™ 96 Well Polycarbonate PCR Microplate, Model M (#6511)	
	Corning® 384 Well Polypropylene PCR Microplate (#6500)	
	Corning® 384 Well Polyethylene PCR Microplate (#6502)	

Order Information

Part #	Description
Thermal Cycler	
7013001	Q-Cycler II Server Standard Thermal Cycler including block – please specify standard block required
7013002	Q-Cycler II Server Gradient Thermal Cycler including block – please specify gradient block required
7013003	Q-Cycler II Server Ultra Gradient Thermal Cycler including block – please specify ultra block required
7013004	Q-Cycler II Satellite Standard Thermal Cycler including block – please specify standard block required
7013005	Q-Cycler II Satellite Gradient Thermal Cycler including block – please specify gradient block required
7013006	Q-Cycler II Satellite Ultra Gradient Thermal Cycler including block – please specify ultra block required
Thermal Blocks	
7004001	Standard Block for 96x0.2ml/96well plate
7004002	Standard Block 96x0.2/48x0.5ml
7004003	Block for 4 Slides/Microarrays
7004004	Gradient Block for 96x0.2ml/96well plate
7004005	Gradient Block 96x0.2/48x0.5ml
7004006	Gradient Block for 384 well plate
7004007	Ultra Gradient Block for 96x0.2ml/96well plate
7004008	Ultra Gradient Block for 384 well plate
Manual	
QBTCMAN	Quanta Biotech Thermal Cycler Manual Operator manual for the thermal cycler and the thermal cycler Software

Sales & Service

Quanta Biotech Ltd

Unit 4 Byfleet Technical Centre

Canada Road

Byfleet

Surrey

KT14 7JX

United Kingdom

Phone +44 (0) 1932 344550 • Fax +44 (0) 1932 353108

Email: info@quantadirect.com

Web: www.Quantabiotech.com

Index

- AC Power Connection, 11
- Adding New User Accounts, 19
- Adding or Modifying User Accounts, 19
- Administrator Login, 18
- Administrator Settings, 15
- Automatic Hot Start command, 23
- Barcode Input command, 23
- Block Exchange, 45
- Changing Users, 32
- Cleaning, 41
- Close Lid command, 24
- Connecting Auto-Q Satellite, QB-96 and QB-24 Thermal Cyclers to the Auto-Q Server, 10
- Connecting Optional Devices, 11
- Creating PCR Programs, 21
- Cycler Management, 15
- Deleting a Program Step, 30
- Deleting Programs, 31
- DND Programming, 21
- Editing Cycler Data, 16
- Editing Programs, 29
- Electrical, 6
- End Cycle command, 26
- First start-up, 18
- Functional Overview, 9
- Fuses, 11
- Gradient Step command, 26
- Gradient Step Conversion, 30
- Grounding Requirements., 12
- Initial Inspection, 8
- Installation, 6
- Loading the Cycler, 34
- Maintenance, 41
- Manual Hot Start command, 24
- Modifying User Accounts, 19
- Networking the Thermal Cycler, 10
- Operating Power Requirements, 11
- Operation, 32
 - Changing Users, 32
 - Opening and Closing the Lid (QB-24 and QB96), 33
 - Running a Program, 35
 - Sample Loading, 34
 - Selecting Programs, 33
 - Setting the Lid Pressure, 35
 - Terminating a Run, 35
 - User Login, 32
 - Viewing Script Temperature Profiles
 - Prior to Run, 33
- Order Information, 51
- Package Contents, 8
- Program Execution, 35
- Program Steps, 23
 - Automatic Hot Start, 23
 - Barcode Input, 23
 - Close Lid, 24
 - End Cycle, 26
 - Gradient Step, 26
 - Manual Hotstart, 24
 - Script Information, 27
 - Start Cycle, 26
 - Store, 26
 - Temperature Step, 24
 - Touchdown, 25
- Program Wizard, 27
- Programming, 21
 - Converting a Gradient Step, 30
 - Deleting a Program Step, 30
 - Deleting Programs, 31
 - Editing Programs, 29
 - General Guidelines, 28
 - Program Wizard, 27
- Reports, 37

- Deleting Reports, 38
- Viewing Reports, 37
- Requirements, 6
- Running a Program
 - Starting a Run, 35
- Safety
 - Instrument ~, 2
 - Symbols and Conventions, 2
- Safety Warnings, 2
- Sales & Service Centers, 52
- Script Information command, 27
- Setting Time and Date, 16
- Site, 6
- Software Operation, 13
 - Required Computer Skills, 13
- Specifications, 46
 - Compatible Barcodes, 49
 - Compatible Plate Formats, 50
 - Networking & Barcode Options, 48
 - Power & Dimensions, 48
 - Reports and Validation, 47
 - Thermal Engine Characteristics, 48
- User & File Management, 47
- User Interfaces, 47
- Start Cycle command, 26
- Start Screen, 13
- Starting a Run, 35
- Store command, 26
- Switching Off the Instrument, 39
- Symbols and Conventions, 2
- Temperature Step command, 24
- Terminating a Run, 35
- Thermal Block, 7
- Touchdown command, 25
- Trouble Shooting
 - If a Power Failure Occurs, 42
- Trouble-Shooting, 42
- UnPacking the Thermal Cycler, 8
- USB Memory Stick Operation, 40
- User Accounts
 - Deleting Accounts, 20
- User Levels, 17
- User Login, 32
- User Management, 17