


# Leica S4 E with A-Stand

Congratulations on the purchase of your new StereoZoom®. Soon you will be enjoying some of the finest imaging capabilities modern technology can provide. All items are packed individually to secure optimal protection for storage and shipping. Prior to unpacking or assembly please make sure that your shipment contains the following individual inner boxes:

## Kit # 10 810 014


- 1 Box 10446293 S4 E StereoZoom
- 1 Box 10447130 10x Eyepiece, fix
- 1 Box 10447131 10x Eyepiece, adj.
- 1 Box 10446339 Focus column
- 1 Box 10446340 Base
- 1 Box 10447039 Dust cover

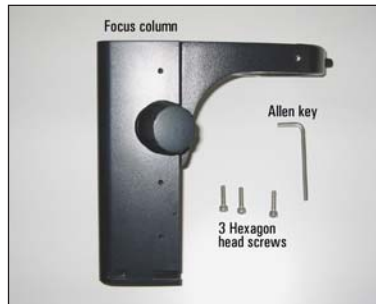

Open your box 10446399 "Focus column" and you will find:




Open box 10446340 "Base"

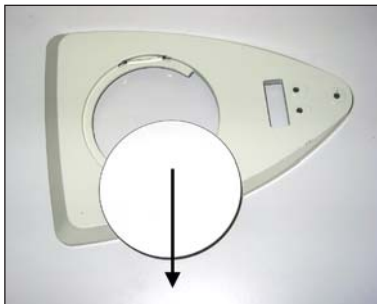

Remove stage plate for assembly of the focus column



Insert 3 screws from the bottom of the base and screw them tightly

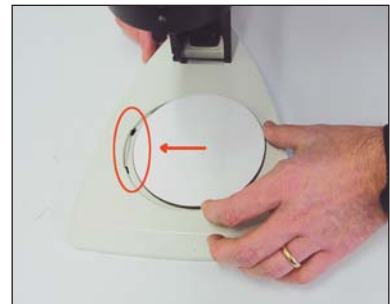

Slide in the stage plate from the side as shown in the picture



Unpack 10446293 StereoZoom® Leica S4 E



Insert your StereoZoom® ...



... in the focus column



Fix StereoZoom® with clamping screw by turning screw clockwise



Remove tube caps from the StereoZoom®



Unpack the eyepiece 10447130 and the eyepiece 10447131



Insert eyepiece 10447130 into the left eyetube as far as it will go



Insert eyepiece 10447131 into the right eyetube as far as it will go



Your StereoZoom® S4 E is now assembled and ready to use

Replacement parts and most common accessories:

Eyeguards	10 447 149
10x eyepiece fix	10 447 130
10x eyepiece adj.	10 447 131
16x eyepiece fix	10 447 132
16x eyepiece adj.	10 447 133
20x eyepiece fix	10 447 134
20x eyepiece adj.	10 447 135
Lens shield	10 446 324

**For care instructions please read your user manual on page 34**

**Please read the user manual for use and safety instructions before putting your StereoZoom® into operation.**

You will enjoy your StereoZoom even more with proper illumination. We have designed the Leica L2 illumination system exclusively for your StereoZoom®. If you want to find out more please visit our web site [www.stereozoom.com](http://www.stereozoom.com) or call us in the US toll free at 800 248 0123.

If you have any questions or problems related to your new StereoZoom®, please call our US toll free number 800 248 0123 and then follow the instructions to reach our customer care service.

**Leica**

MICROSYSTEMS