

# Installation and Operating Instructions for the FluidX IntelliCode software

Rev 1, release date: July 2014



FluidX Ltd. Monks Heath Hall, Chelford Road, Nether Alderley, Cheshire, SK10 4SY, UK.  
T: +44(0) 1625 861 614 F: +44(0) 1625 861 615 E: [info@fluidx.eu](mailto:info@fluidx.eu) [www.fluidx.eu](http://www.fluidx.eu)

This document is for information only; the manufacturer accepts no liability for errors contained herein or for incidental or consequential damages with the furnishing, performance, or use of this material.

Unless otherwise specified references to names or parts is purely casual and has the purpose of illustrating the product. The contents of this publication may not be reproduced in any form or by any means (including electronic storage and retrieval or translation into a foreign language) without prior agreement and written consent from the copyright owner.

The information contained in this document is subject to change without notice.

#### **Table of Contents:**

- 1. Welcome**
- 2. Introduction**
- 3. Installation**
- 4. Activation**
- 5. Deactivation**
- 6. Managing licenses**
- 7. Getting Started**
- 8. Exporting Files**
  - Text**
  - Document**
  - Destination and Header Information**
  - Database**
  - Audit**
- 9. Selecting an instrument**
- 10. Creating a new profile**
  - No reads**
  - Selecting FluidX 2d Rack ID**
  - Selecting SBS Rack Barcode**
- 11. Features**
  - Toolbar**
  - Virtual Imager**
  - Export Options**
  - Rack and tube format**
  - Orientation**
  - Rack id**
  - SBS Sample rack**
- 12. Decode Process**
  - Parameters**
  - Crop**
  - Image Process**
  - Validate**
  - No tube**
  - Decoder**
- 13. Remote Command Set**

## **Welcome:**

?????

## **Introduction:**




Intellicode™ is designed to have no or very little user input. Profiles have been pre-installed for the majority of consumable types.


## **Installation:**

**Do not connect a reader device until installation is complete.**

Ensure the user installing the software has full admin rights, (Read, Write and Modify, to the registry)


Close all other applications during the installation process

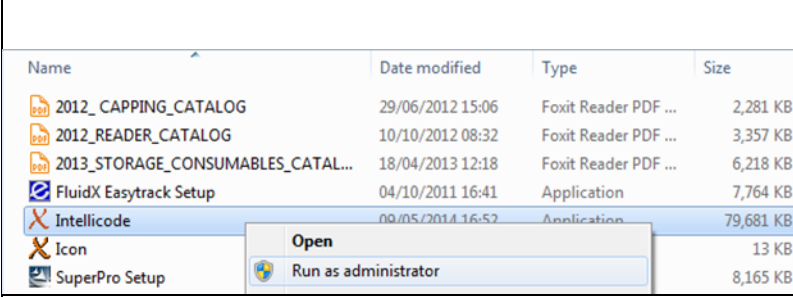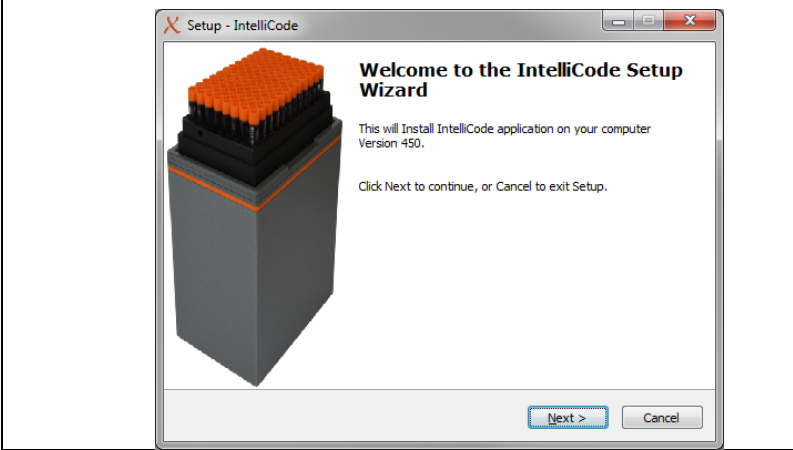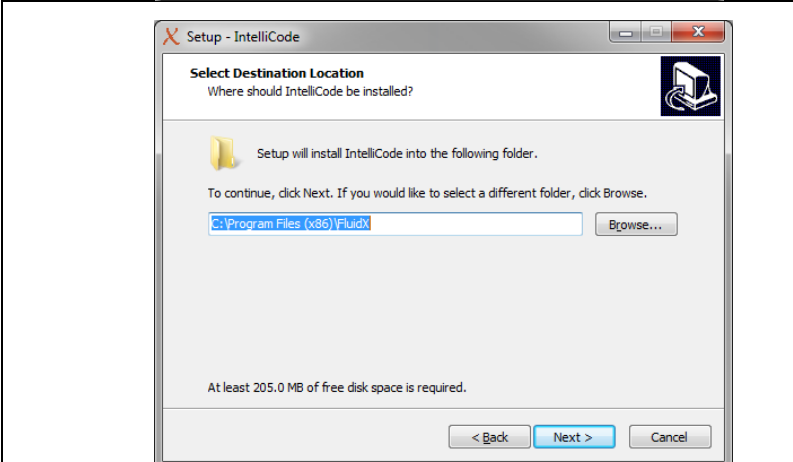	<p>Open the IntelliCode folder on the fluidx software USB stick.</p> <p>Run the IntelliCode application as administrator</p>
	<p>Follow the on screen instructions.</p> <p>Select "Next" at the Welcome screen to start the install process.</p>
	<p>Verify the Install path and click "Next"</p> <p>The install will proceed.</p>

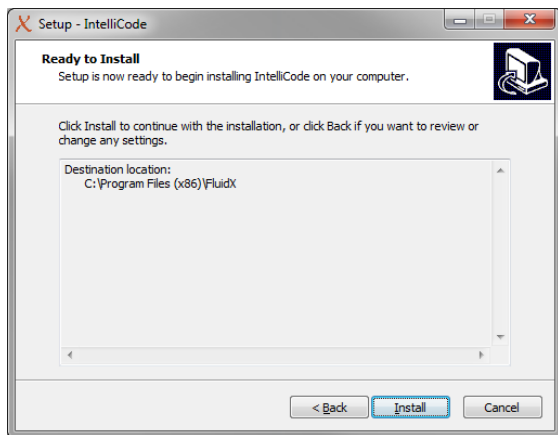


Verify the Install Details and click "Install"

The install will proceed

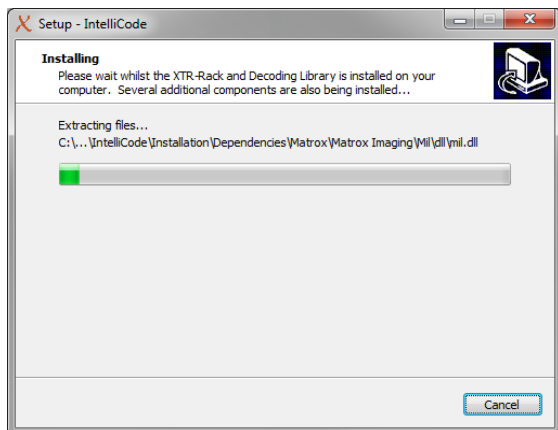


The installation will proceed through various steps of the installation.

This may take a few moments

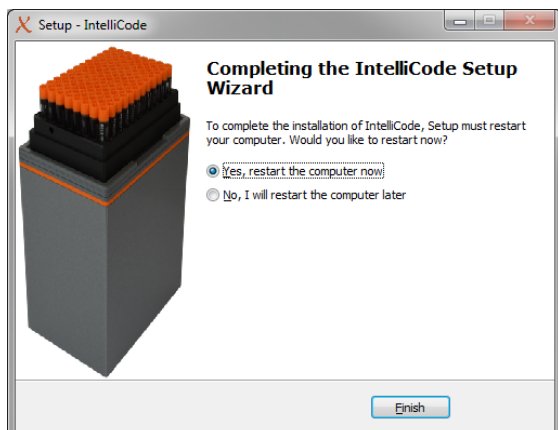

Select "Yes, restart the computer now", then select "Finish"

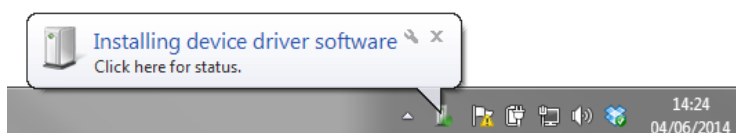


Once the restart has been performed, plug the scanning device into an available port on the PC.

Wait for the device to be recognised.




IntelliCode has now been successfully installed on the PC, to open the software double click the shortcut on the desktop.

## Activation:

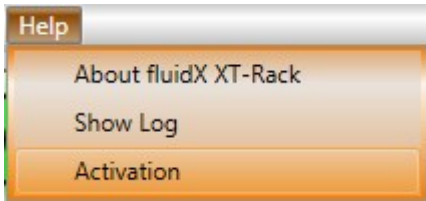


From the Tool bar select "Help" then "Activation"

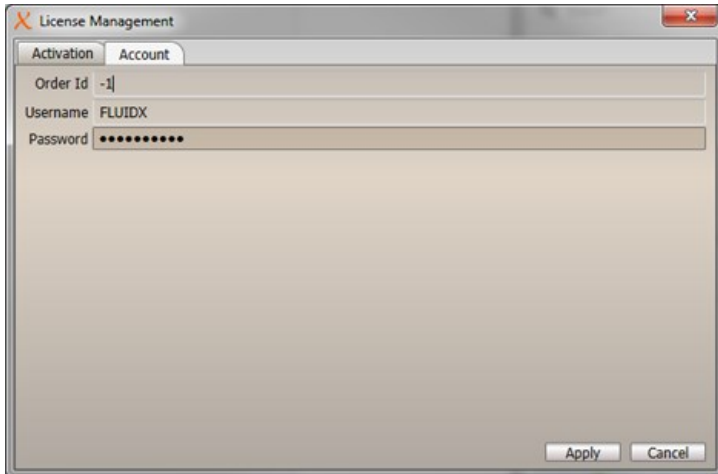

Go to the "Account" tab and enter your Username and Password.

Select Apply

Account information, order Id, Username and Password will be included with the software package.  
(Credit card/ note pad on install disk, information on USB install sticker)

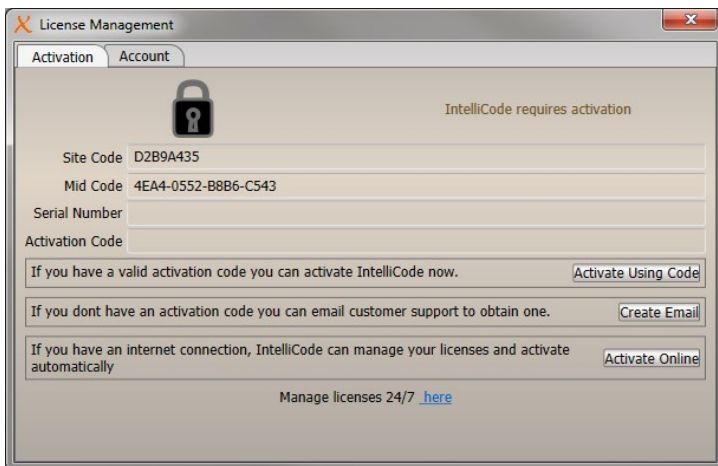


The software can be used for 7 days without activation, however after 7 days the software will become locked.

Activation can be performed at any time after installation.

To activate your IntelliCode software go to "Help" and select "Activation"


The options below show how to activate the software.

1. Activate on Line, internet connection needed.

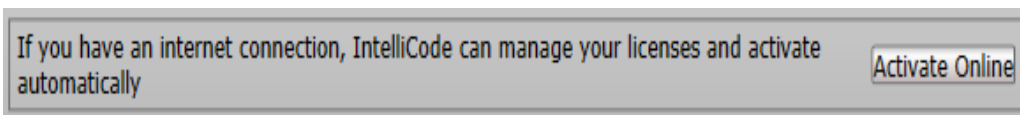

Activation is complete.

## 2. Create Email

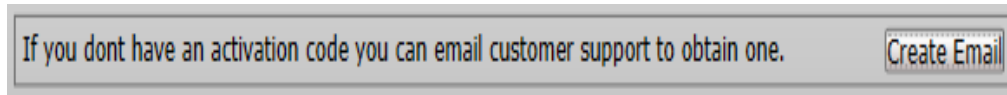


Automatically creates an email with the necessary information to send to the FluidX customer support. Replies should be received within 24 hours. Once the activation code is received from customer support, follow step three below.

## 3. Activate using code

If the serial number and activation number is already known, enter these into the fields on the activate screen.

Serial Number	DBF3-45A2-BF-5D0E-DEMO
Activation Code	0C8DFADF-B823D6B9-118E8165-B9B2DEMO

Select "Activate using code"



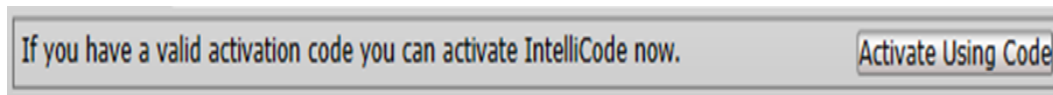


Activation is now complete.

## **Deactivation:**

Unless additional licenses are purchased the software can only be activated on one PC at a time. If the license needs to be migrated to a different PC, this can be achieved by deactivating and transferring the license.

This can be managed via the Activation screen.

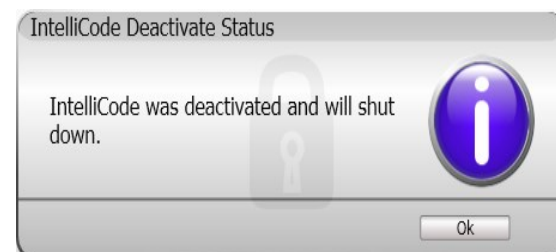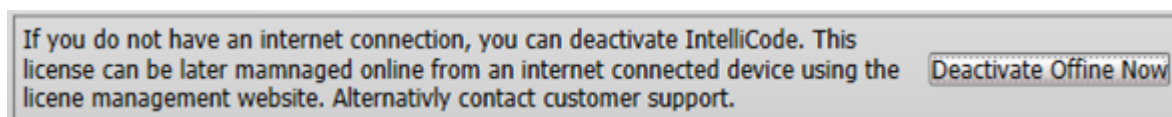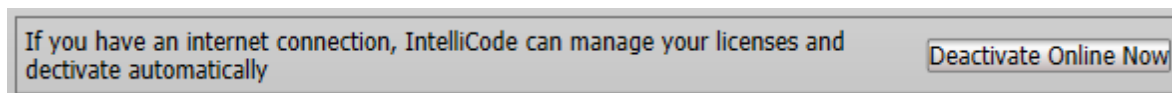

If the software is needed in a new location, then once the install has taken place, follow the activation steps in section ???

## **Managing Licenses**

Licenses can be controlled via the on-line activation centre, this may be more useful in the instance when there are additional licenses assigned to one scanner, or multiple scanners are installed on site.

Access can be gained by selecting the "Manage licenses 24/7" or going to [www.fluidxactivation.co.uk](http://www.fluidxactivation.co.uk)

Manage licenses 24/7 [here](#)

## fluidX activation centre

Username:

Password:

☒ Remember me

Login

[Forgot password?](#)

Enter the Username and Password.

This can be found on **credit card** received with the software or on text note on the installation USB.

If the PC is not shared and in a safe location "Remember me" can be selected.

Select the order number by pressing the key icon.

An additional box will appear below with activation information regarding that order.

For a new activation select the red Plus sign and enter the Site and mid code from the IntelliCode activation screen.

The Note field can be used for internal use, for example which computer holds the licence, or department.

## fluidX activation centre


Logged as: FLUIDX








Account

Orders

Log out










### ORDERS AND LICENSES

This is the list of your active orders. Please click on  button to view licenses for selected order.

	ID	Program name	Date	Licenses	Payment type
	81800	Fluidx Xtr-flex pro	12/12/2013	10 / 3	
	81802	Xtr96	16/12/2013	10 / 0	
	81803	Xtr96	16/12/2013	20 / 1	
	81804	Xtr96	16/12/2013	8 / 0	
	81806	Xtr Single	17/12/2013	3 / 2	
	81822	XtRack	05/02/2014	10 / 4	
	81843	Xtr Single	28/04/2014	5 / 1	

1

Status: 4 of 10 licenses currently activated for this order.

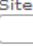	Site code	Machine ID	Note	Activated	Expires	
	D474458C	17DD-102F-EE7B-9E40	xt:Rob	12/03/2014	N/A	
	E614582D	6056-0E34-ED0F-22D9	xt:FLUIDX	13/03/2014	N/A	
	5D8DB2AD	5569-C8DE-06B5-3004	xt:Matt	28/03/2014	N/A	
	36C70845	EF7A-5740-8030-8F31	xt:Rob	16/05/2014	N/A	
						

Site code

D2B9A435

Machine ID (MID)

4EA4-0552-B8B6-C543

Note

Sophie

Add license

Status: Please enter Site code and Machine ID values for new license.

### Select Add licence

Selecting the Email icon option generates a pre-populated email which can be sent to the user, IT department, manager etc. as a record of activation.

The activation code can also be located in the body of text.



This can be used for "Activate using code" if the other activation methods are not available.

**SEND EMAIL (CLIENT: FLUIDX)**

Subject:

Recipient:

CC:

Body:

Program information:  
-----  
Program title: **XtRack**

Computer information:  
-----  
Site code: **D289A435**  
MID code: **4EA4-0552-B886-C543**

License information:  
-----  
Activation date: 02/06/2014  
Validity: Unlimited  
Note: Sophie  
Activation code:  
**188287E1-A2FD13A5-2C4FEA32-1893C359**  
-----  
fluidX Support.


**Send**

Status: **Email successfully sent.**

## Getting Started:

IntelliCode is designed to have no or very little user configuration.  
Profiles have been pre-installed for the majority of consumable types.

Open the IntelliCode software by using the shortcut on the desktop, or from the IntelliCode folder.  
C:\Program Files (x86)\FluidX\IntelliCode

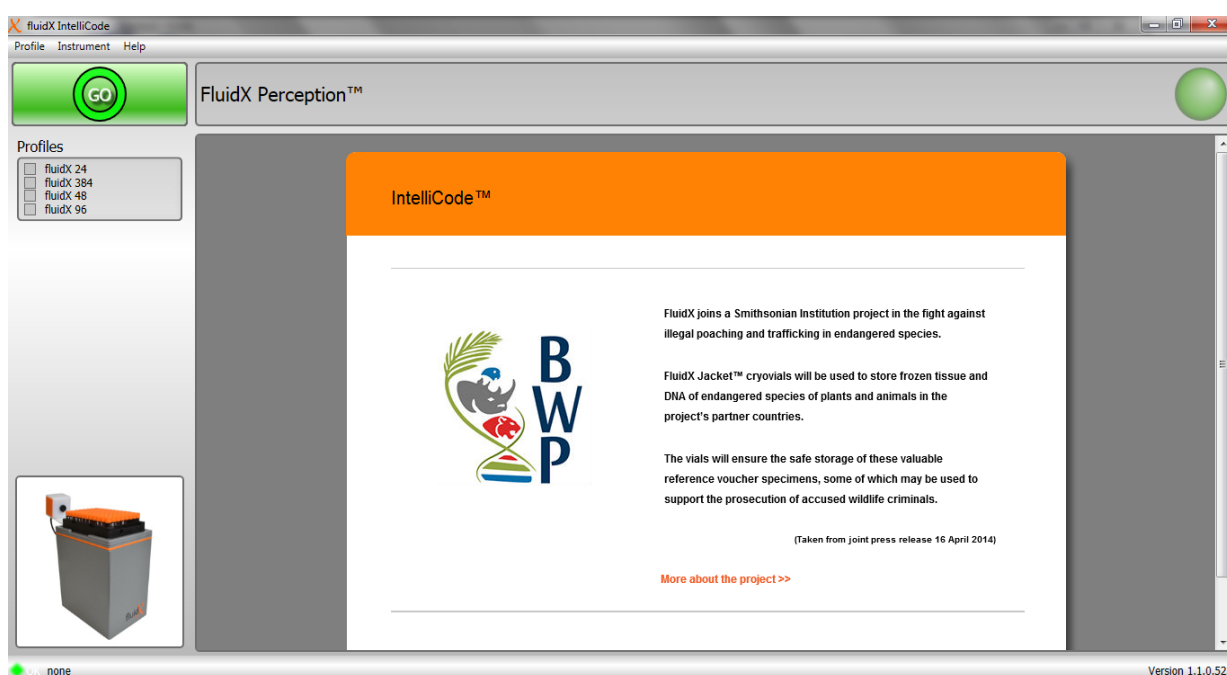

Select an in-

**strument.** See section 9: Selecting an Instrument


Place a rack on the reader and select which profile matches the manufacturer and configuration of the rack. If the manufacturer or configuration is not known, select all the profiles.

Select "GO"




The grid box positions will turn green when all the tubes are decoded.



The number of successful, failed and no Tube output is report at the top of the grid with decode time.

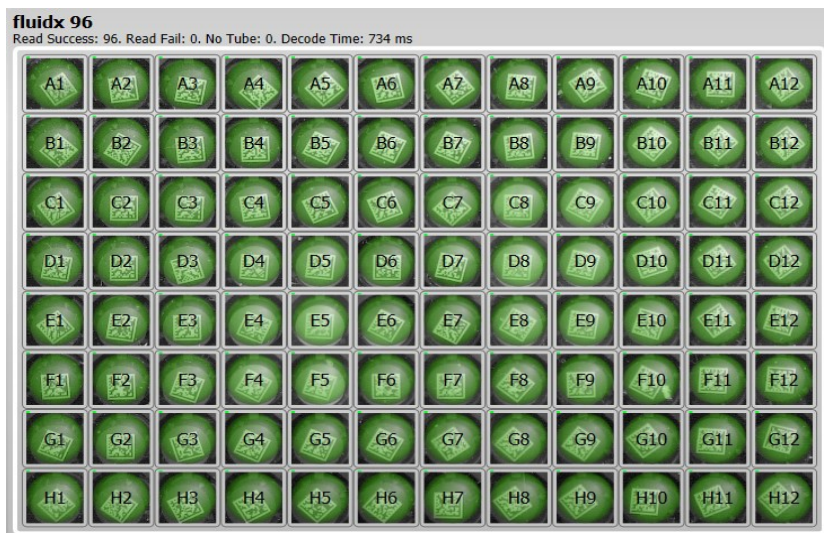

Click on a single cell to see more detailed information of the code.

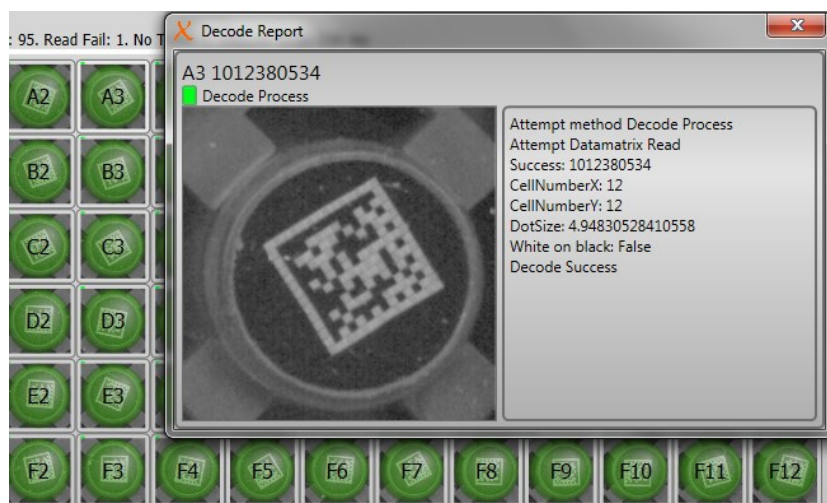

Click on the 1d\_rackid\_1 for more detailed information of the linear barcode.

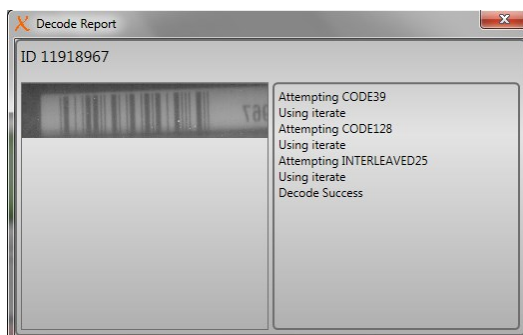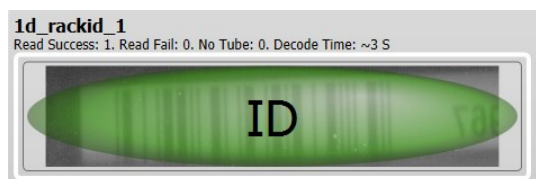


Below the grid table is the destination directory for the exported data.

Right clicking on the Export line gives the options

**Open Containing Folder:** Opens the Export destination folder

**Open With Associated Application:** Open the Export data file.

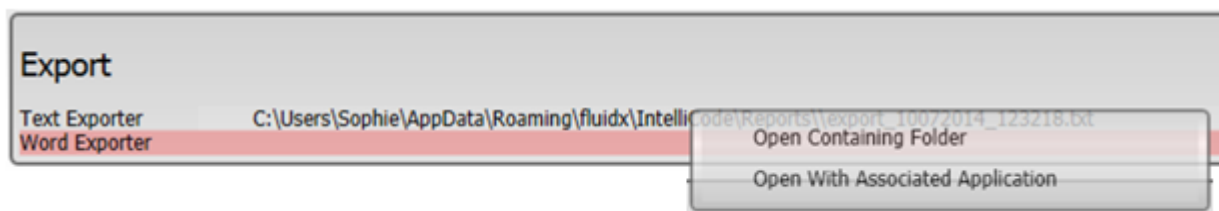


Session Success/Fails are below the Export data path.

Session Success/Fails				
Success				Fail
fluidx 96 .A1 FC12798817	fluidx 96 .A2 FC12798825	fluidx 96 .A3 FC12798833	fluidx 96 .A4 FC12798841	
fluidx 96 .A5 FC12798849	fluidx 96 .A6 FC12798857	fluidx 96 .A7 FC12798865	fluidx 96 .A8 FC12798873	
fluidx 96 .A9 FC12798881	fluidx 96 .A10 FC12798889	fluidx 96 .A11 FC12798897	fluidx 96 .A12 FC12798905	
fluidx 96 .B1 FC12798818	fluidx 96 .B2 FC12798826	fluidx 96 .B3 FC12798834	fluidx 96 .B4 FC12798842	
fluidx 96 .B5 FC12798850	fluidx 96 .B6 FC12798858	fluidx 96 .B7 FC12798866	fluidx 96 .B8 FC12798874	
fluidx 96 .B9 FC12798882	fluidx 96 .B10 FC12798890	fluidx 96 .B11 FC12798898	fluidx 96 .B12 FC12798906	
fluidx 96 .C1 FC12798819	fluidx 96 .C2 FC12798827	fluidx 96 .C3 FC12798835	fluidx 96 .C4 FC12798843	
fluidx 96 .C5 FC12798851	fluidx 96 .C6 FC12798859	fluidx 96 .C7 FC12798867	fluidx 96 .C8 FC12798875	
fluidx 96 .C9 FC12798883	fluidx 96 .C10 FC12798891	fluidx 96 .C11 FC12798899	fluidx 96 .C12 FC12798907	

In the event a tube cannot be read, go to section 10: Creating a new profile, No reads




## Exporting Files

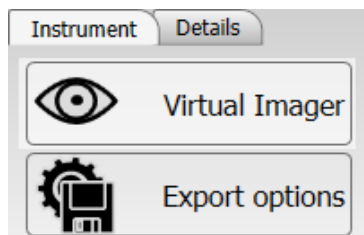

To access the Export options, double click the profile.

To configure your exported files, select Export options

### Text:

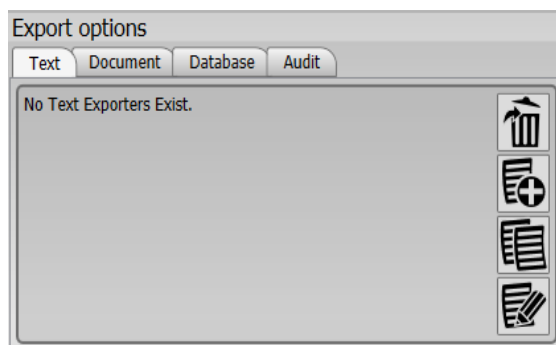

Delete



Delete an Export type

Create a new Export


Copies the Export

Edit the Export

The export file can be given a custom name by typing in the space by "Name"

Auto Create automatically imports the evaluators.


The export information can be customised by copying and pasting the selection from Evaluators to the Edit column. Changing the extension to .csv creates an Excel file

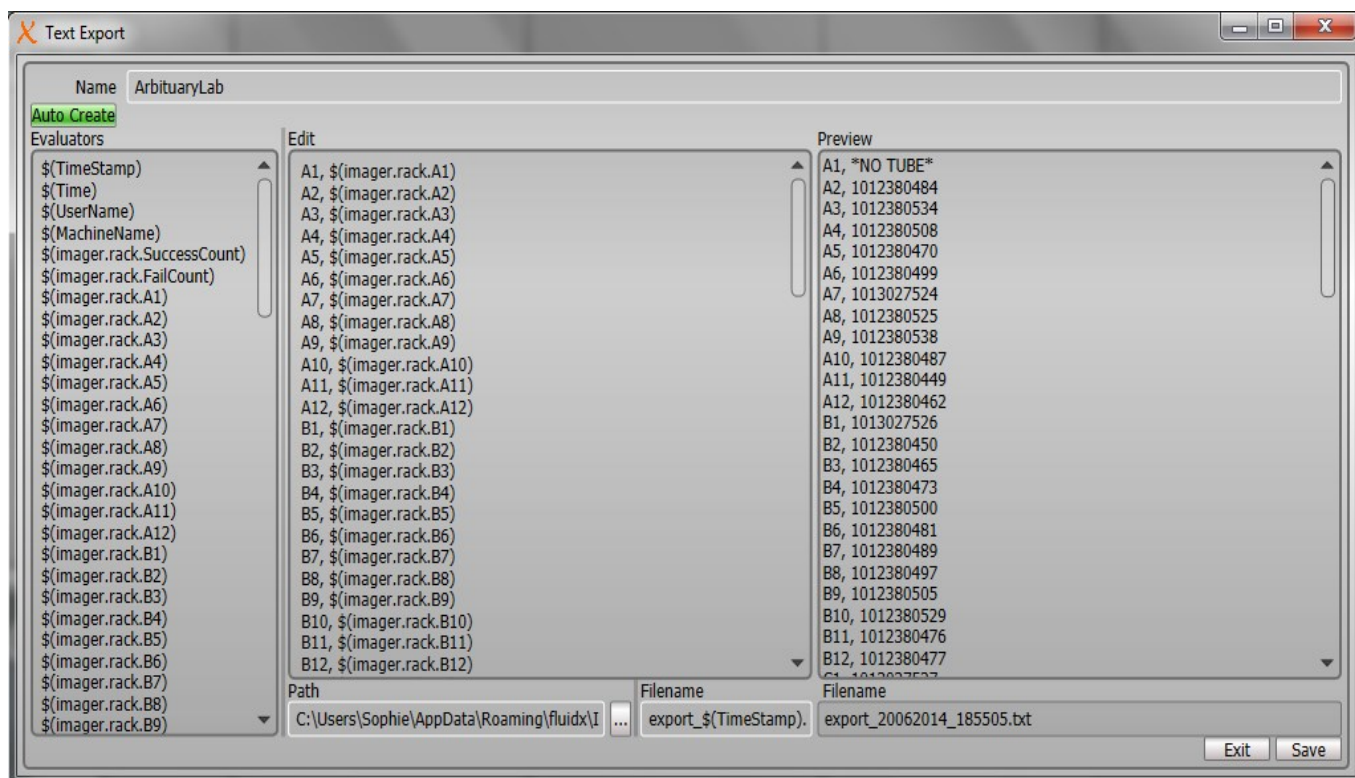




## Document:

To create a document file select Document from the Export Option and click to Create a new Word Exporter

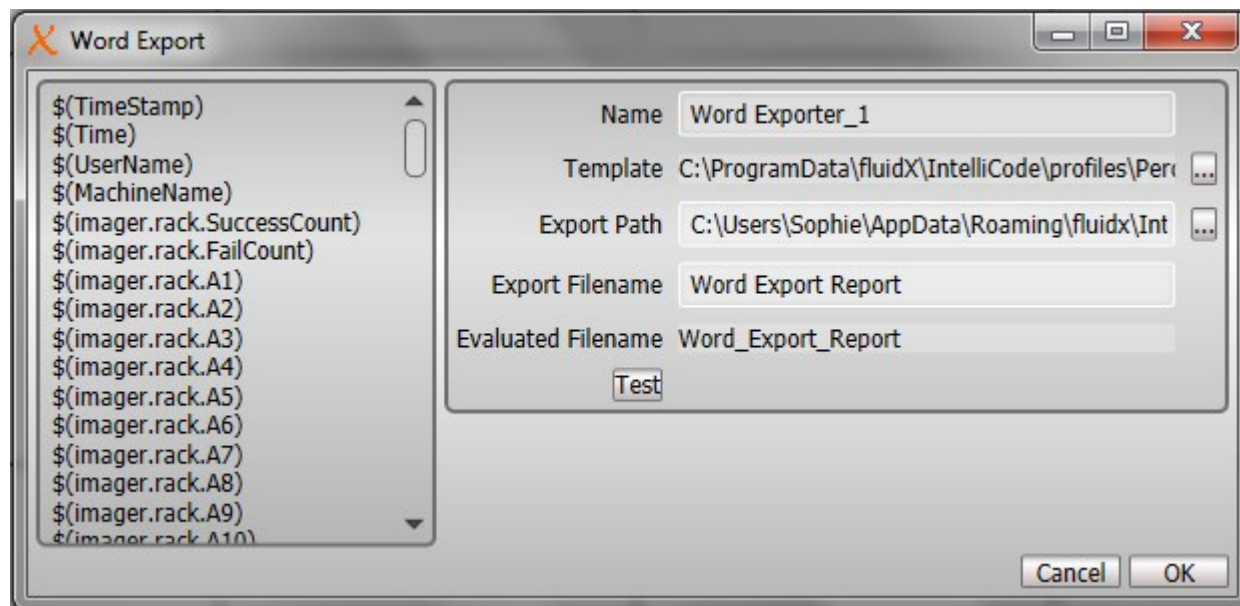

The 'Word Export' dialog box contains a list of variables on the left and configuration fields on the right. The variables list includes: \$(TimeStamp), \$(Time), \$(UserName), \$(MachineName), \$(imager.rack.SuccessCount), \$(imager.rack.FailCount), \$(imager.rack.A1) through \$(imager.rack.A10). The configuration fields on the right are: Name (Word Exporter\_1), Template (C:\ProgramData\fluidX\IntelliCode\profiles\Perception), Export Path (C:\Users\Sophie\AppData\Roaming\fluidX\IntelliCode), Export Filename (Word Export Report), and Evaluated Filename (Word\_Export\_Report). A 'Test' button is located below the Evaluated Filename field. At the bottom right are 'Cancel' and 'OK' buttons.

Templates have been pre-configured and stored at C:\ProgramData\fluidX\IntelliCode\profiles\Perception. Select "Test" for an example of the output, select "OK" to save and migrate to the Export page.

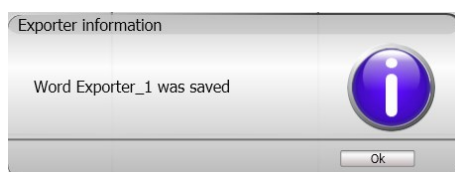

## IntelliCode Word Export Example

### Standard 96 Report Template


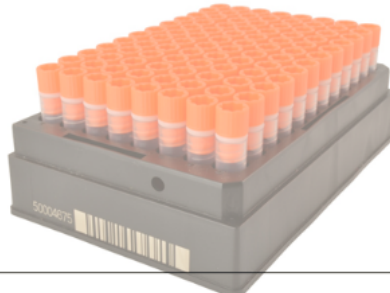
ComputerID: MachineName

UserID: UserName









Number of Tubes Decode Fails: decodefails

Rack Id: RackId

	1	2	3	4	5	6	7	8	9	10	11	12
A	A1	A2	A3	A4	A5	A6	A7	A8	A9	A10	A11	A12
B	B1	B2	B3	B4	B5	B6	B7	B8	B9	B10	B11	B12
C	C1	C2	C3	C4	C5	C6	C7	C8	C9	C10	C11	C12
D	D1	D2	D3	D4	D5	D6	D7	D8	D9	D10	D11	D12
E	E1	E2	E3	E4	E5	E6	E7	E8	E9	E10	E11	E12
F	F1	F2	F3	F4	F5	F6	F7	F8	F9	F10	F11	F12
G	G1	G2	G3	G4	G5	G6	G7	G8	G9	G10	G11	G12
H	H1	H2	H3	H4	H5	H6	H7	H8	H9	H10	H11	H12





## Example Header Information:

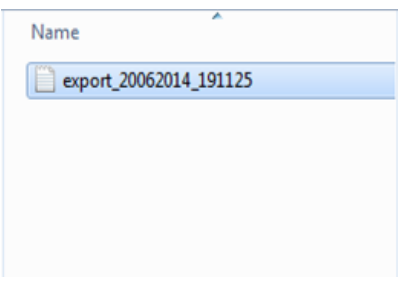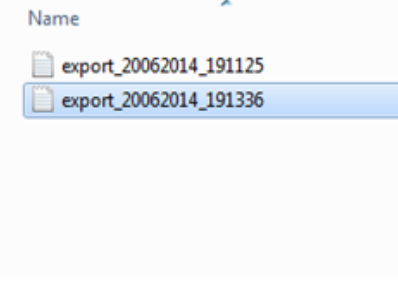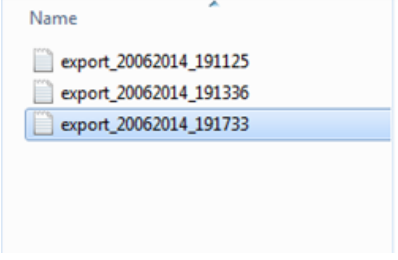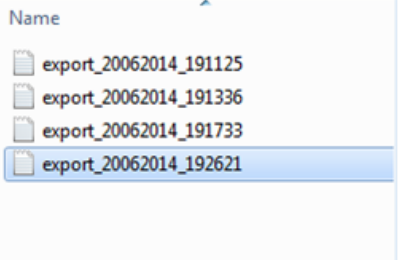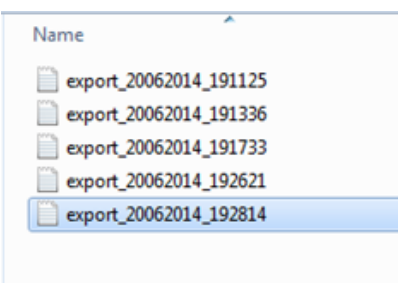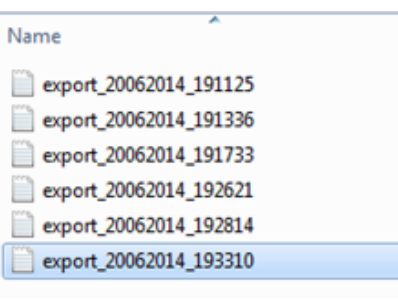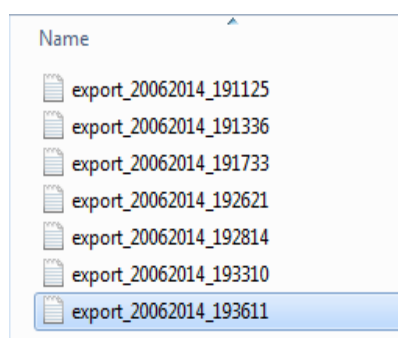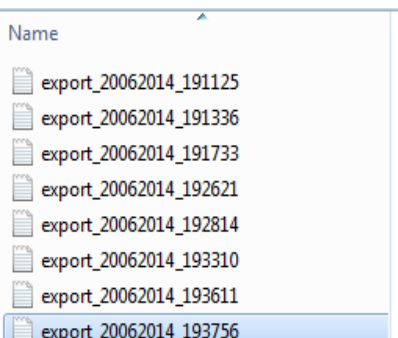No Header information	Time Stamp (Date_Time)
 <p>             Name              A1, *NO TUBE*              A2, 1012380484              A3, 1012380534              A4, 1012380508              A5, 1012380470              A6, 1012380499              A7, 1013027524              A8, 1012380525              A9, 1012380538              A10, 1012380487              A11, 1012380449              A12, 1012380462           </p>	 <p>             Name              20062014_191336              A1, *NO TUBE*              A2, 1012380484              A3, 1012380534              A4, 1012380508              A5, 1012380470              A6, 1012380499              A7, 1013027524              A8, 1012380525              A9, 1012380538              A10, 1012380487              A11, 1012380449              A12, 1012380462           </p>
Time	User Name
 <p>             Name              20/06/2014 19:17:33              A1, *NO TUBE*              A2, 1012380484              A3, 1012380534              A4, 1012380508              A5, 1012380470              A6, 1012380499              A7, 1013027524              A8, 1012380525              A9, 1012380538              A10, 1012380487              A11, 1012380449              A12, 1012380462           </p>	 <p>             Name              Sophie              A1, *NO TUBE*              A2, 1012380484              A3, 1012380534              A4, 1012380508              A5, 1012380470              A6, 1012380499              A7, 1013027524              A8, 1012380525              A9, 1012380538              A10, 1012380487              A11, 1012380449              A12, 1012380462           </p>
Machine Name	Successful decode number (imager.rack.SuccessCount)
 <p>             Name              SOPHIE-THINK              A1, *NO TUBE*              A2, 1012380484              A3, 1012380534              A4, 1012380508              A5, 1012380470              A6, 1012380499              A7, 1013027524              A8, 1012380525              A9, 1012380538              A10, 1012380487              A11, 1012380449              A12, 1012380462           </p>	 <p>             Name              95              A1, *NO TUBE*              A2, 1012380484              A3, 1012380534              A4, 1012380508              A5, 1012380470              A6, 1012380499              A7, 1013027524              A8, 1012380525              A9, 1012380538              A10, 1012380487              A11, 1012380449              A12, 1012380462           </p>
Failed decodes (imager.rack.FailCount)	All selected
 <p>             Name              0              A1, *NO TUBE*              A2, 1012380484              A3, 1012380534              A4, 1012380508              A5, 1012380470              A6, 1012380499              A7, 1013027524              A8, 1012380525              A9, 1012380538              A10, 1012380487              A11, 1012380449              A12, 1012380462              B1, 1013027526           </p>	 <p>             Name              20062014_193756              20/06/2014 19:37:56              Sophie              SOPHIE-THINK              95              0              A1, *NO TUBE*              A2, 1012380484              A3, 1012380534              A4, 1012380508              A5, 1012380470              A6, 1012380499              A7, 1013027524              A8, 1012380525              A9, 1012380538              A10, 1012380487           </p>

## Legacy Rack Enabled:

Exports the data in the same configuration as the rack.

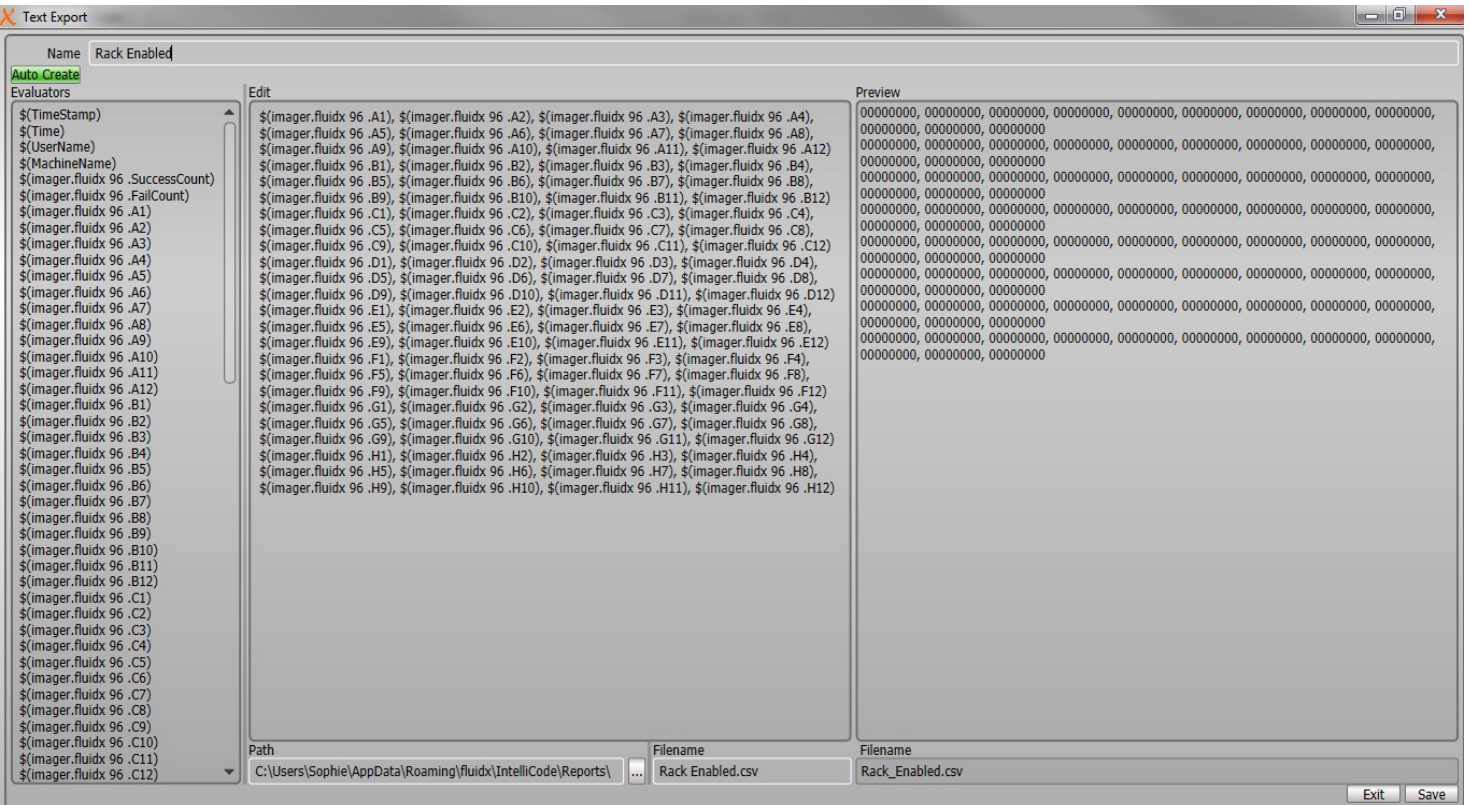

Enter a comma for the value to be in the next column.

Enter a return for the data to be in the row below.

P10

✕

✓

fx

	A	B	C	D	E	F	G	H	I	J	K	L	M
1	FC12798817	FC12798825	FC12798833	FC12798841	FC12798849	FC12798857	FC12798865	FC12798873	FC12798881	FC12798889	FC12798897	FC12798905	
2	FC12798818	FC12798826	FC12798834	FC12798842	FC12798850	FC12798858	FC12798866	FC12798874	FC12798882	FC12798890	FC12798898	FC12798906	
3	FC12798819	FC12798827	FC12798835	FC12798843	FC12798851	FC12798859	FC12798867	FC12798875	FC12798883	FC12798891	FC12798899	FC12798907	
4	FC12798820	FC12798828	FC12798836	FC12798844	FC12798852	FC12798860	FC12798868	FC12798876	FC12798884	FC12798892	FC12798900	FC12798908	
5	FC12798821	FC12798829	FC12798837	FC12798845	FC12798853	FC12798861	FC12798869	FC12798877	FC12798885	FC12798893	FC12798901	FC12798909	
6	FC12798822	FC12798830	FC12798838	FC12798846	FC12798854	FC12798862	FC12798870	FC12798878	FC12798886	FC12798894	FC12798902	FC12798910	
7	FC12798823	FC12798831	FC12798839	FC12798847	FC12798855	FC12798863	FC12798871	FC12798879	FC12798887	FC12798895	FC12798903	FC12798911	
8	FC12798824	FC12798832	FC12798840	FC12798848	FC12798856	FC12798864	FC12798872	FC12798880	FC12798888	FC12798896	FC12798904	FC12798912	
9													
10													
11													
12													


## Destination and header information:

**Path:** Destination path for the data, selecting the “....” option allows the user to browse the directory.  
*Default path* = %appdata%\fluidx\IntelliCode\Reports





Quick access to the folder can be achieved by clicking on the "FluidX Data" shortcut on the desk top.

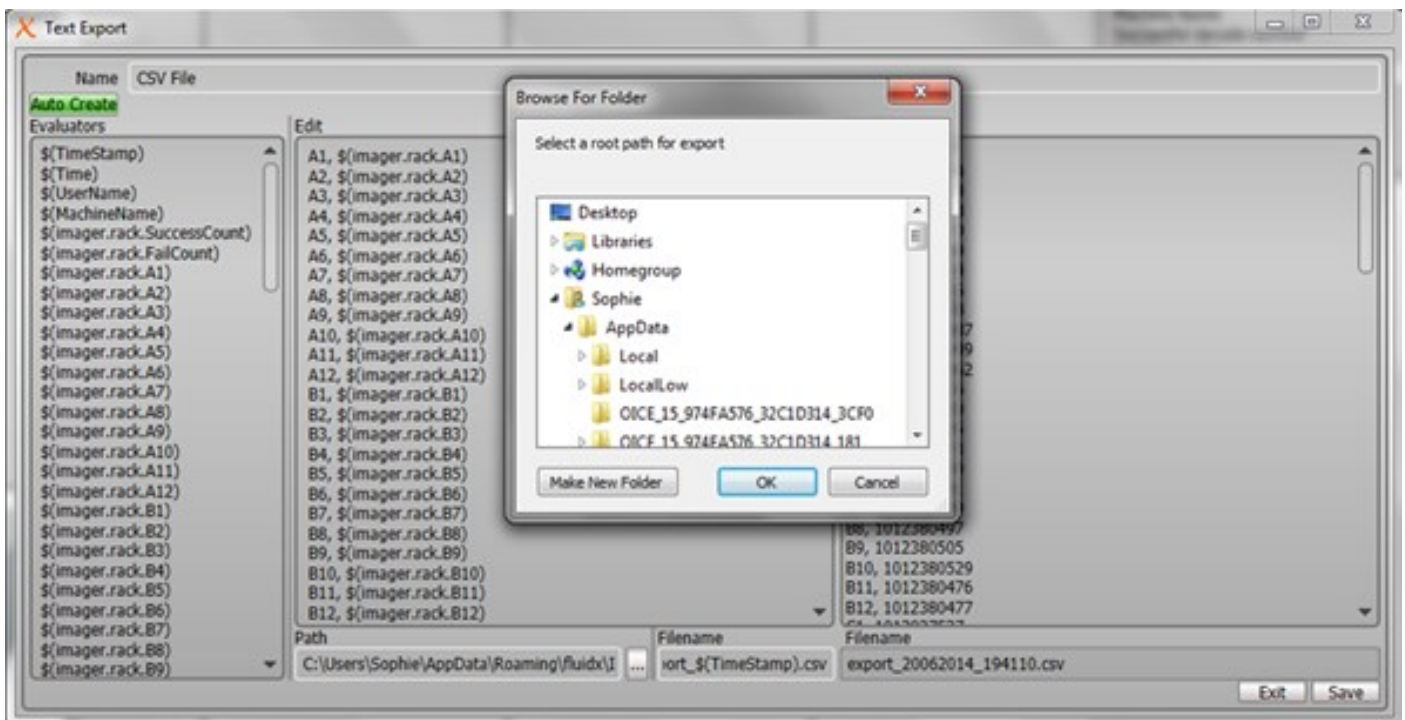


**Filename:** Sets the name of the file when saved


**Filename:** Example of how the file name will look when exported.

## Database


Database allows data transfer with ODBC

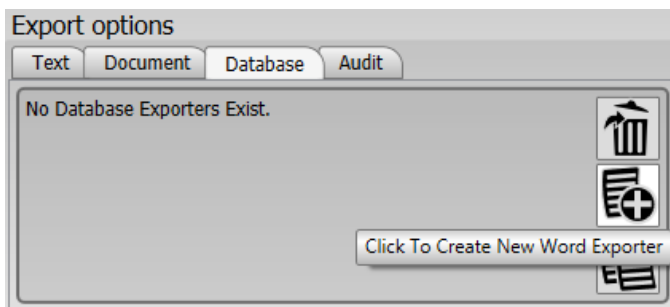

From the Export options section select "database" and select New Word Exporter

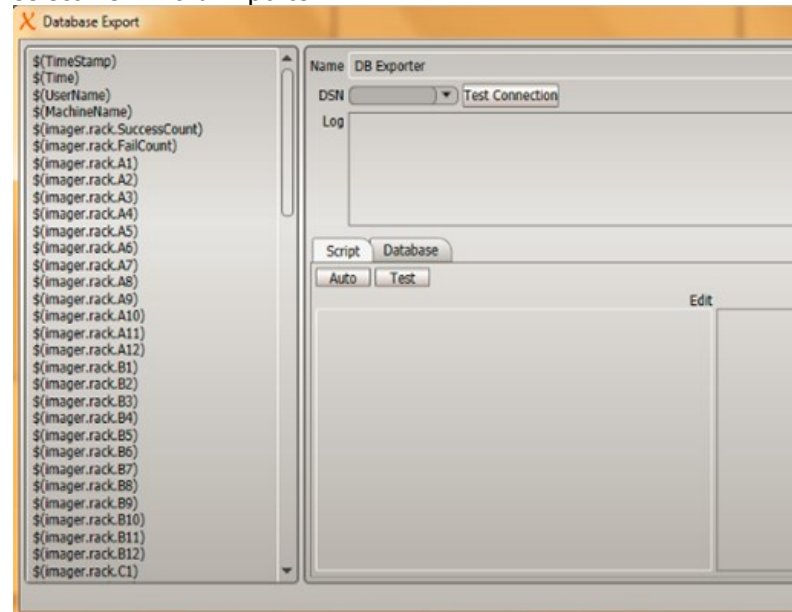



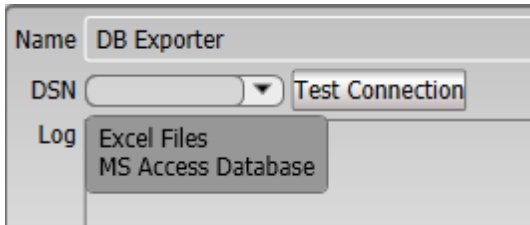


From DSN select the compatible log.

Using either copy and paste or "Auto" move the required information to the Edit column.  
An example of the export is displayed in the Preview column.  
Select "Save"




### Audit

The Audit option allows scanned files to be retrieved if the Export destination has not been previously created.

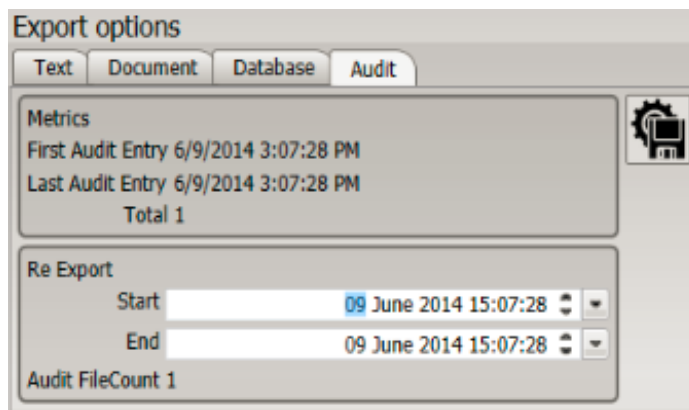

## Selecting an Instrument

IntelliCode is compatible with all fluidx readers.

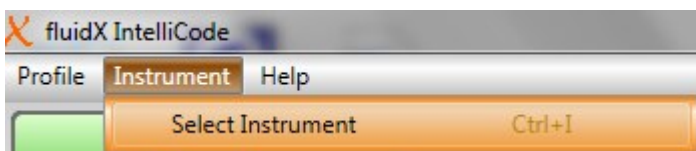


Go to Instrument and "Select Instrument"

Various readers can be chosen from a list of instruments, if the text is greyed out, the device has not been installed, or not installed correctly.


### Virtual instrument:

This option can be used in the instance when an integrator or prospect purchaser would like to try the software

Select "Show All", click on reader type and press "Select"  
Making a selection will show to the left a description of the device.

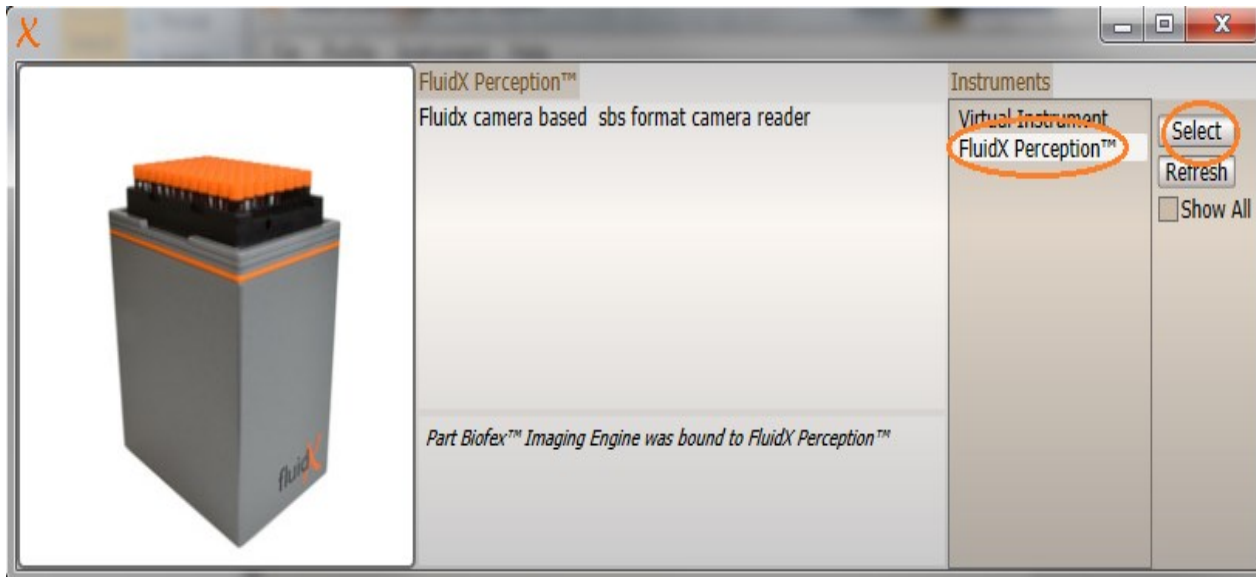


## Creating a new profile

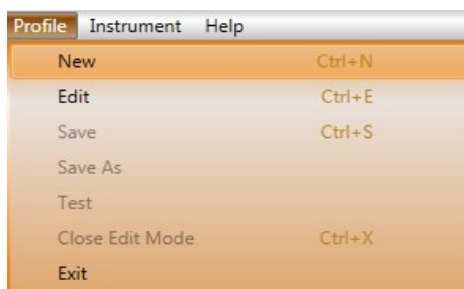


Go to "Profile" and Select "New"

Using the mouse drag the desired "SBS Sample Rack" over the rack image

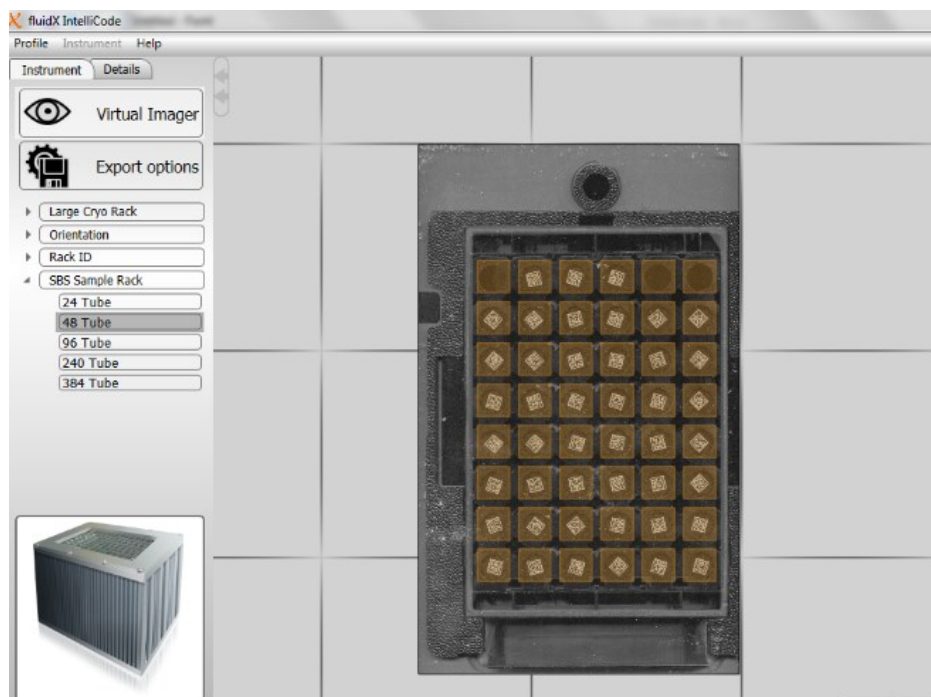

The aim is to match the centre quadrant rather than the whole rack.

Refine the Region of Interest (ROI). This improves decode time and optimises decoding for difficult to read codes.

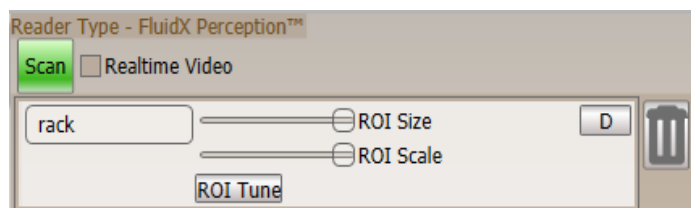

"rack"- Ability to change the rack name.



"ROI Tune" Automatically centres and optimises the ROI

"D" Returns the setting to factory default.

Select "Decode"




If there are no tubes present then it will be necessary to set up the no tube/read feature. *See Section 12 Decode Process; No Tube.*

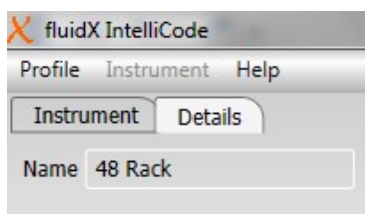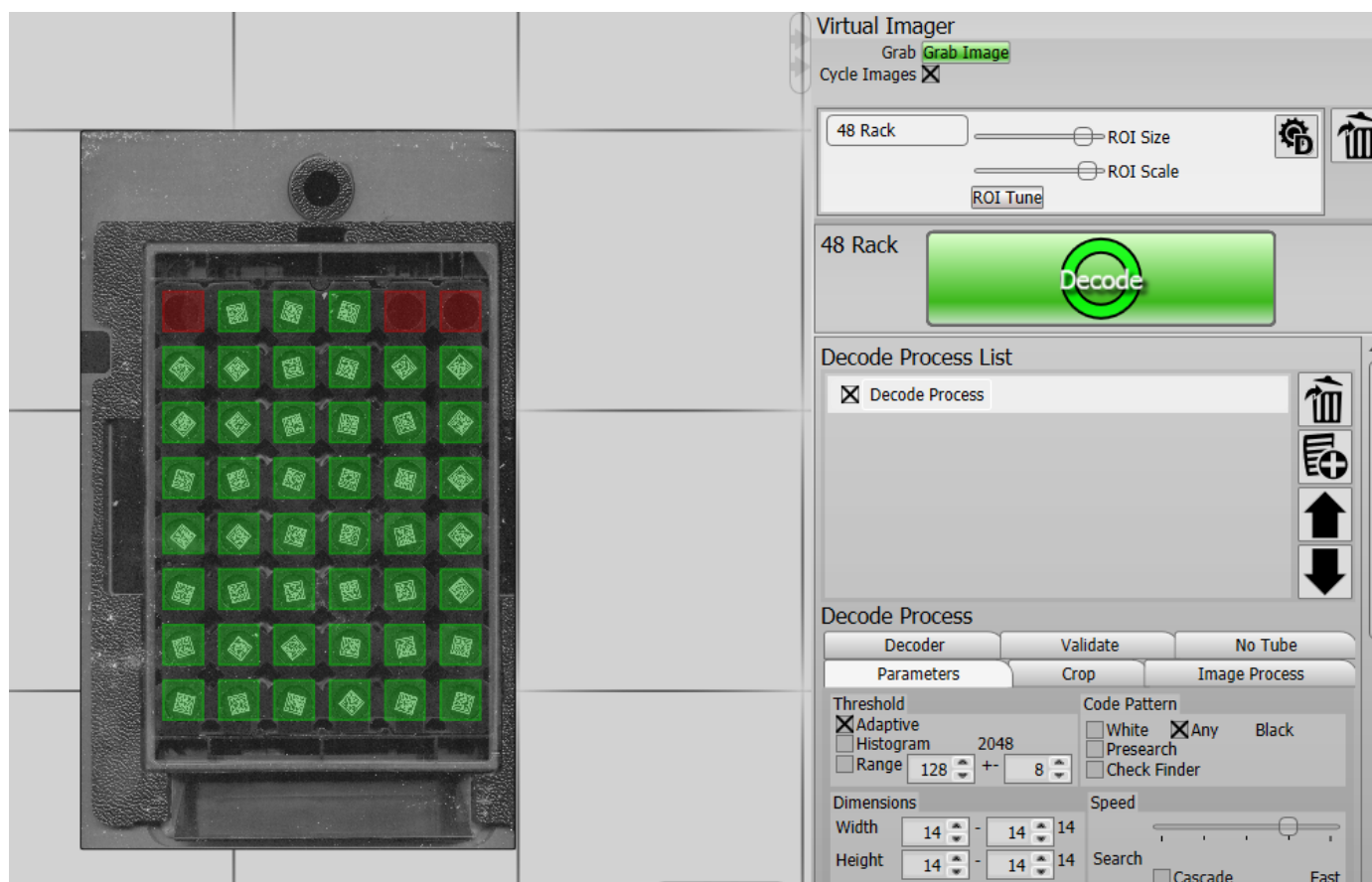


The Name of the rack can be generated and saved by going to the Details tab


To save the new profile go to "Profile" the "Save As"

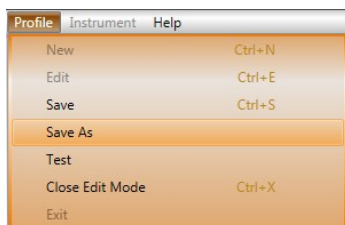

Default folder: C:\ProgramData\fluidX\IntelliCode\profiles\Virtual Instrument

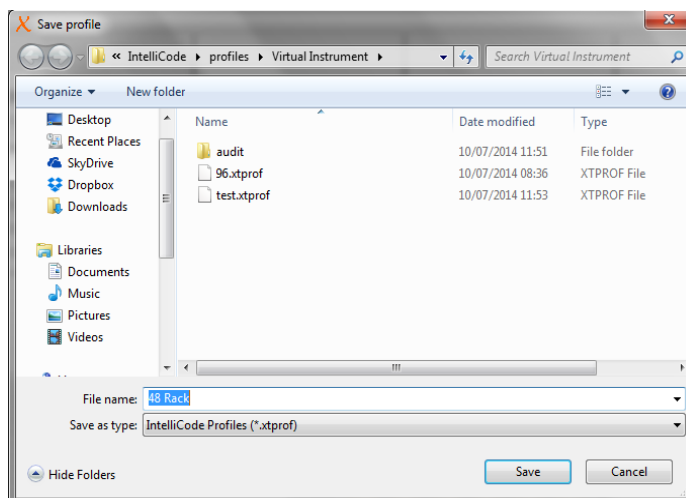


Once the file is saved, go to "Profile" then "Close Edit Mode".

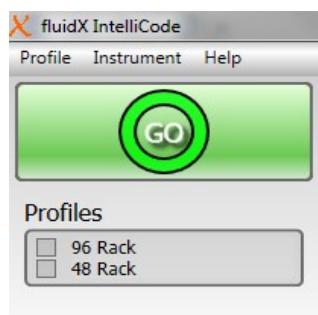

The new profile will be added to the default list on the main screen.


To select the profile, check the desired box.

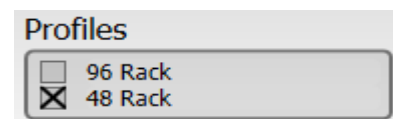

Multiple profiles can be selected and IntelliCode will automatically detect the rack configuration.

## Selecting FluidX 2d Rack ID


Expand Rack ID and drag FluidX 2d Rack ID over the image of the ID tube.

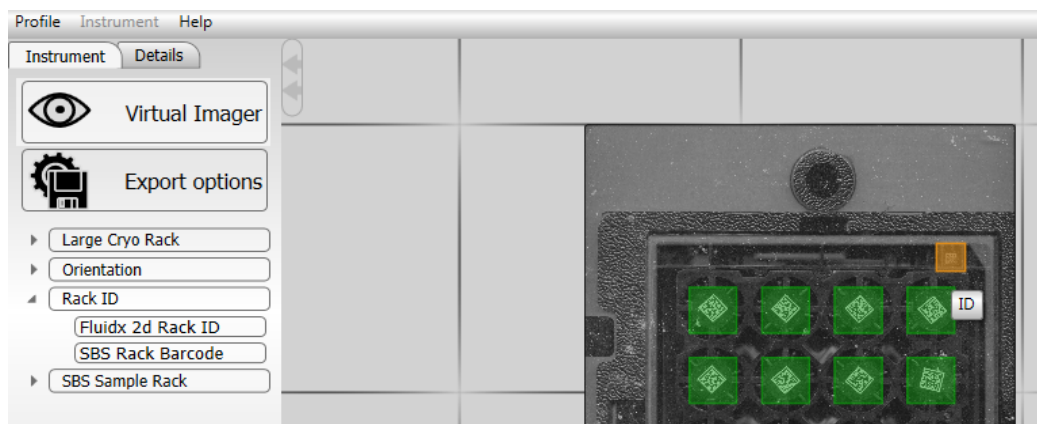

Select Decode




Scroll down in the Decode Process list to see the image of the ID code.

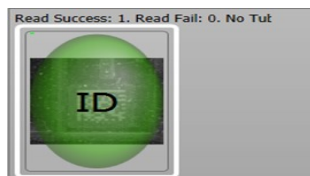

Clicking on the icon generates the decode report




In the event that the 2d\_rack\_1.ID cannot be read, a dialogue box will appear so the value can be manually entered.

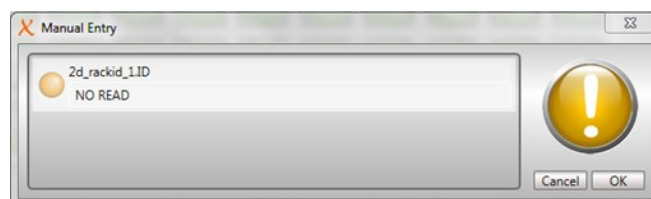

## Selecting SBS


Drag the option for SBS Rack Barcode over the linear barcode.

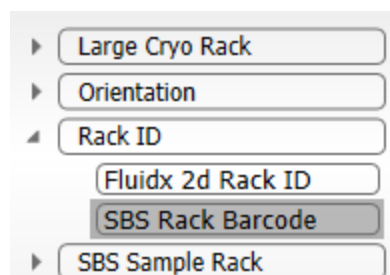

## Rack Barcode



Select 1d\_rackid\_

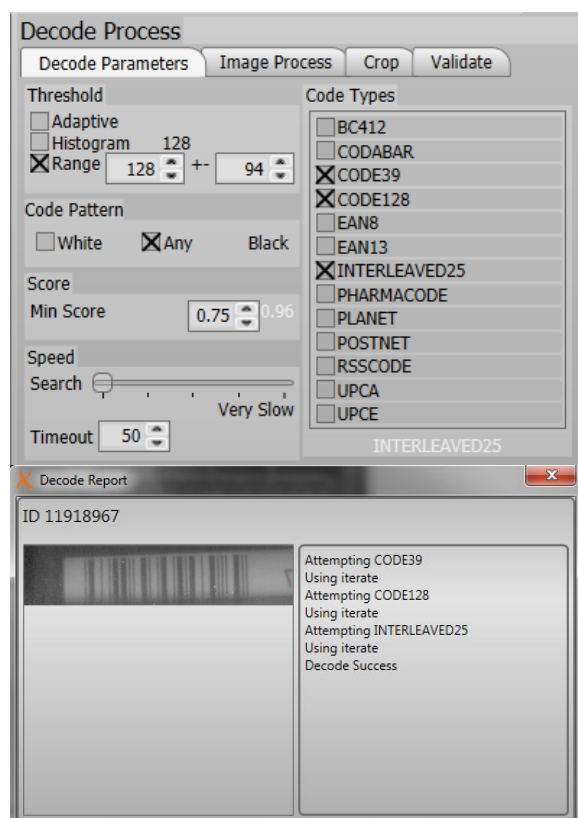

The majority of linear barcodes will CODE39, CODE128 or INTERLEAVED25

Adjustment of the Decode Parameters may be necessary for non-standard codes.

Clicking on the decode image shows a dialogue box with a detailed Decode Report.




## No Reads

If tubes are not being read, the decode process should be checked.

To determine the parameters of a code click on a green cell in the grid under the Decode Process List and compare the parameters to a code that has not been read.


Cell Number, Width and Height and Foreground colour are the most common things to change.  
Check the code is within the ROI and is clear, not damaged and free from frost or debris on the glass.

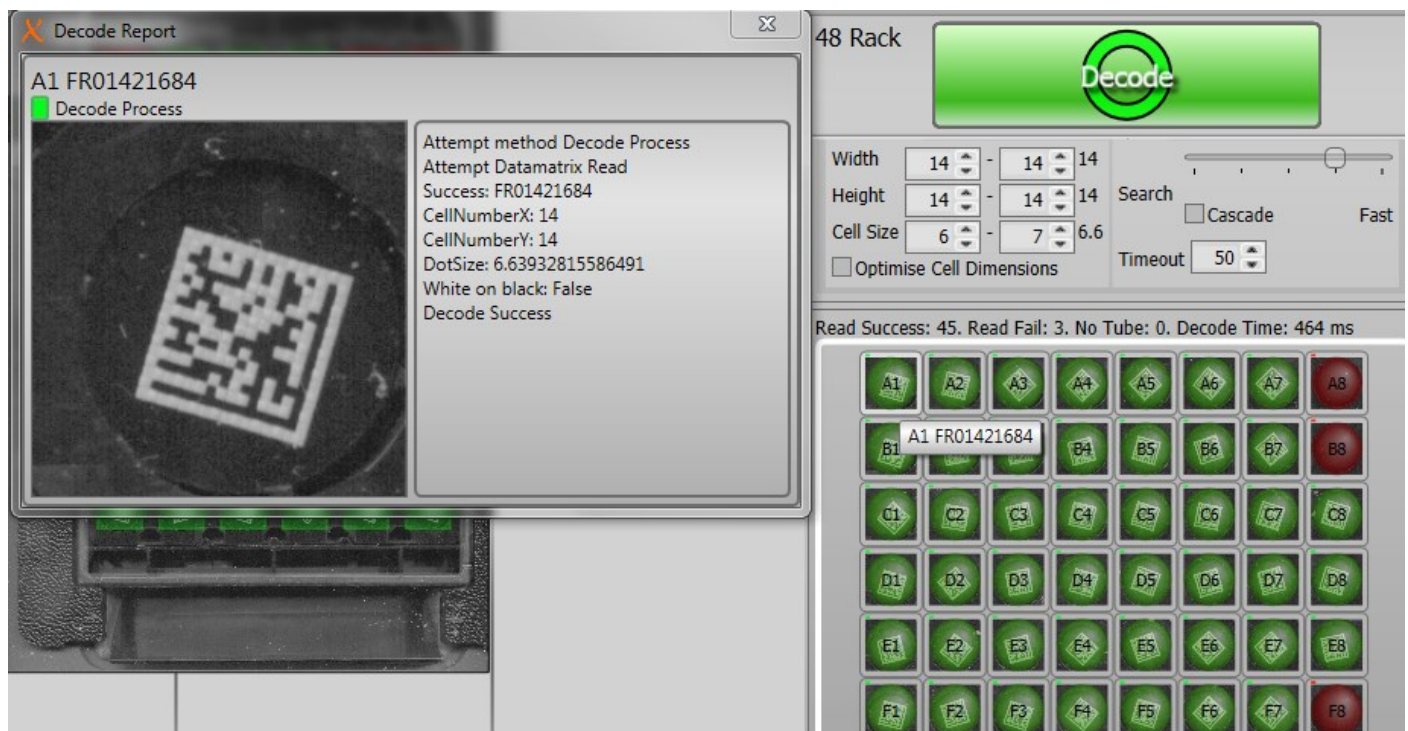

## Features

### Tool bar

#### Profile

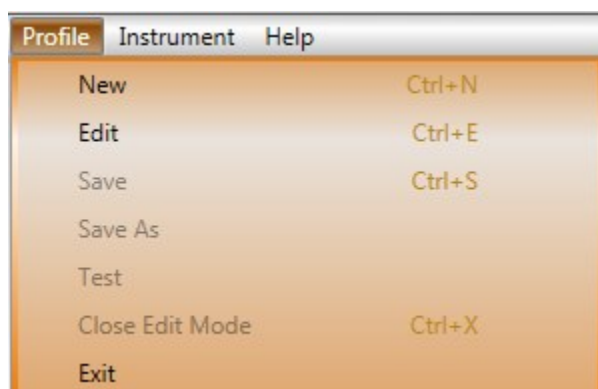

**New** – Creates new rack configuration

**Edit**- Changes can be made to a current configuration

**Save**- saves a xtpo file


**Save As**- saves a xtpo file

**Test**- Performs a decode so the configuration can be checked before closing the edit mode.

**Close Edit Mode**- Saves and closes settings and returns to main page

**Exit-** Closes the software

## Help

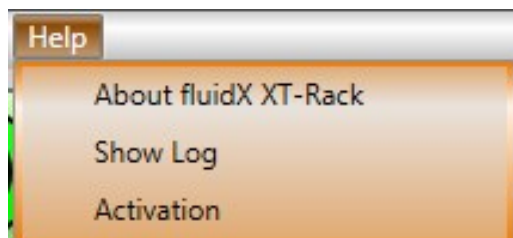


**About fluidX XT-Rack-**Licensing, manufacturer and software information


**Show Log:** Transcript of software activity.

For problems encountered with IntelliCode this should be copied and pasted, and sent to [Support@fluidx.eu](mailto:Support@fluidx.eu)

**Activation:** Information on IntelliCode activate and account settings. See section: **Activation**

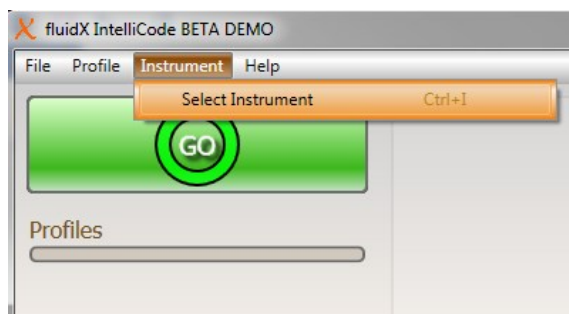


**Instrument:** Shows installed devices

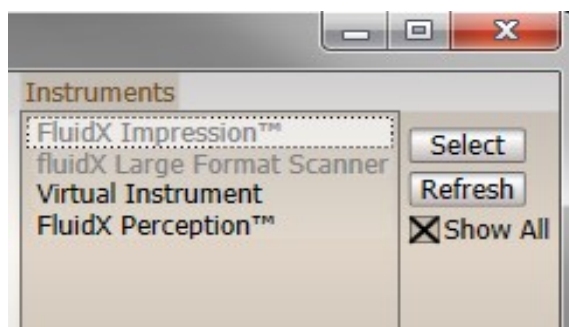



Various readers can be chosen from a list of instruments, if the text is greyed out, the device has not been installed.

Select "Show All", click on reader type and press "Select"  
Making a selection will show to the left a description of the device.

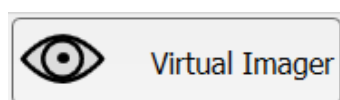


**Virtual Imager:** Allows IntelliCode to work without any hardware. Software can be used for evaluation with a virtual instrument.

A box will appear to the right of the screen.  
Select "Grab Image" scroll through the images until the correct format is shown the screen

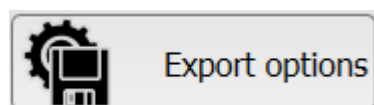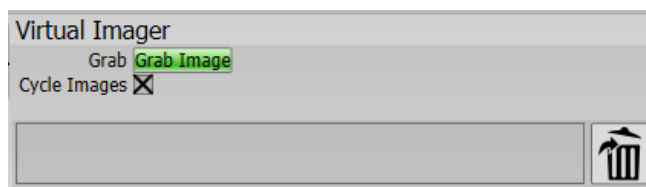

See section 8, Exporting files.

**Export Options:** Allows the user to export the reported values of the codes to various formats such as text file and excel.

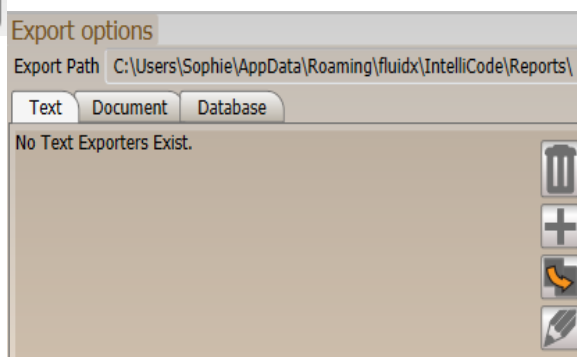

A vertical menu with four options: Large Cryo Rack, Orientation, Rack ID, and SBS Sample Rack. Each option is in a rounded rectangular button.

**Rack and Tube Format:**

A sub-menu for 'Large Cryo Rack' with two options: 81 Tube and 100 Tube. Each option is in a rounded rectangular button.

**Large Cryo Rack:** For use with the FluidX Perception large format reader

A sub-menu for 'Orientation' with two options: 1d orientation and 2d orientation. Each option is in a rounded rectangular button.

**Orientation:** Used to auto-orient the rack.

A sub-menu for 'Rack ID' with two options: Fluidx 2d Rack ID and SBS Rack Barcode. Each option is in a rounded rectangular button.

**Rack ID:** A selection can be made if the Fluidx 2d rack is used or the 1d linear rack ID to record the rack number.

A sub-menu for 'SBS Sample Rack' with five options: 24 Tube, 48 Tube, 96 Tube, 240 Tube, and 384 Tube. Each option is in a rounded rectangular button.

**SBS Sample Rack:** Different rack formats can be selected.

**Decode**  
**Parameters**

The 'Decode Process' dialog box has several tabs: Decoder, Validate, No Tube, Parameters, Crop, and Image Process. The 'Parameters' tab is active. It contains sections for Threshold (Adaptive checked, Histogram unchecked, Range 128 +/- 8), Code Pattern (White unchecked, Any checked, Black unchecked, Presearch unchecked, Check Finder unchecked), Dimensions (Width 12-16-14, Height 12-14-14, Cell Size 3-6-4.1, Optimise Cell Dimensions unchecked), Speed (a slider), Search (Cascade unchecked, Fast selected), and Timeout (50).

**Process**



**Threshold**

Creates a binary image for the decoder. There are a number of methods which achieve this.

**Adaptive:**

Used in situations where the lighting across the image may be uneven and/or the image to decode occupies smaller region of the ROI.

**Histogram:** Used for scanner based readers with a pure black and white image.

**Range:**

Uses a brute force method. Can cause a delay in decoding

**Dimensions**

Information regarding the code to find. The set value to the right is the last recorded pixel value.

**Width:**

The size range for the code width (in dots)

**Height**

The size range for the code height (in dots)

**Cell Size**

The size range for a code dot (in pixels )

The value to the right is a successful value found by the decoder.

The minimum and maximum values should be set to be integers above and below the set value.

**Optimise Cell Dimensions:**

With this option unchecked, the decode process uses the same dimensions for each decode attempt. Checking this option will lock down the decoder to use the dimensions of the first successful decode. If a rack is guaranteed to contain codes of the same dimensions then checking this option can cause a decrease in overall decode time.

Note for racks with mixed code dimension, leave this option unchecked.

**Code Pattern**

Information regarding the code matrix pattern.

**White**

Check this option if the code is white on black. Uncheck for black on white.

**Any**

Check this option if the code can be either black on white or white on black. For racks containing mixed code patterns, ensure this option is checked.

**Presearch**

This should be used when there is a small code within a large ROI area

**Check Finder:** A validation step to prevent code within code detection.


**Speed:** Reducing the speed uses more aggressive algorithms when deciphering a code. Decreasing the speed makes the decoding more robust at the cost of time.

**Cascade:** When checked, each decode attempt will start using the "Decode Method" selected. Retries with successively slower methods will occur until a successful decode has been achieved.

**Time out:** The maximum time ( in milliseconds ) allowed for each decode attempt.

## Crop

Cropping extracts the image which will be used for the decode attempt.

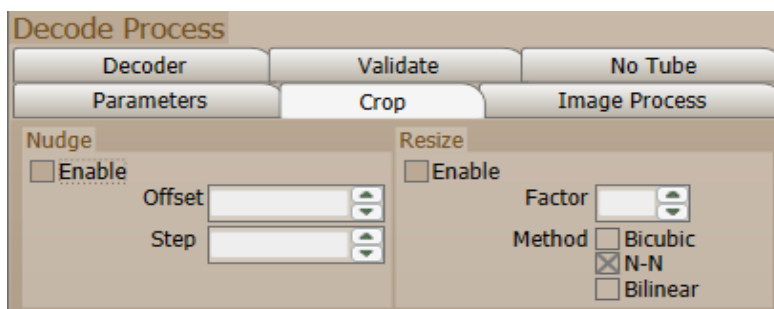

**Nudge:** Offsets the extraction region by a number of pixels. Use when code positions can vary or if codes can fall slightly outside of the ROI.

**Offset:** +/- Maximum offset of region extraction.

**Step:** Iterates from -Offset to +Offset in these steps. Smaller values can increase decode time.

**Re-size:** scales the cropped image. If codes are very small this can improve decode success.

**Factor:** resizes the image by this factor.


**Method:** The scaling method.

Bicubic ( slowest) - [http://en.wikipedia.org/wiki/Bicubic\\_interpolation](http://en.wikipedia.org/wiki/Bicubic_interpolation)

N-N ( fastest) - [http://en.wikipedia.org/wiki/Nearest-neighbor\\_interpolation](http://en.wikipedia.org/wiki/Nearest-neighbor_interpolation)

Bilinear [http://en.wikipedia.org/wiki/Bilinear\\_interpolation](http://en.wikipedia.org/wiki/Bilinear_interpolation)

## Image Process

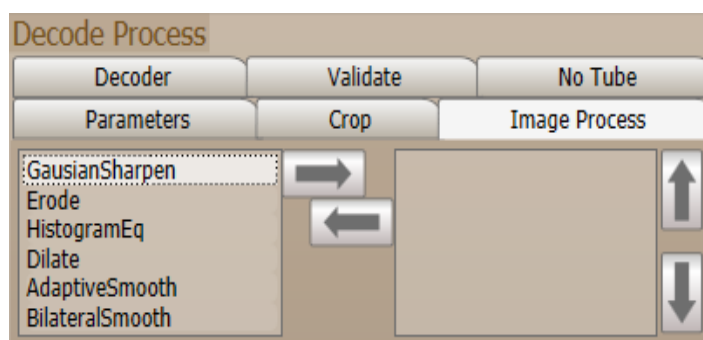

Pre-process the image before attempting decode. This can increase the decode success for damaged, badly marked or lower quality codes.

Image processing works as a stack. Each method is applied to the image from top to bottom.

An overview of the image processing methods are.

**-GaussianSharpen:** Sharpens the image. Can improved decode success with over-marked codes

**-Histogram:** Creates a higher contrast image. Can improved decode success with codes with have little contrast.

**-Dilate:** Attempts to increase the code marking. Can improved decode success with damaged or poorly marked codes.


**-Erode:** Attempts to decrease the code marking. Can improved decode success with over-marked codes.

**-AdaptiveSmooth**

**-BilateralSmooth**

Both smoothing algorithms. Attempts to remove image noise.

## Validate

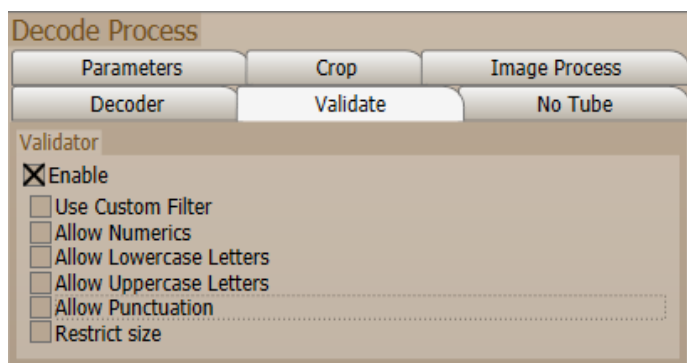


Enabling the Validator expands various code output options.

This stops codes being read inside codes, and allows only codes with letters or numbers

## No Tube

A series of sample lines at varying angles are taken across the cell image and checked for compliance with a set of rules defining a likely matrix code. If the criteria are met the analysis finishes and a matrix code is deemed to be present.

The samples lines can be increased or decreased as deemed necessary.

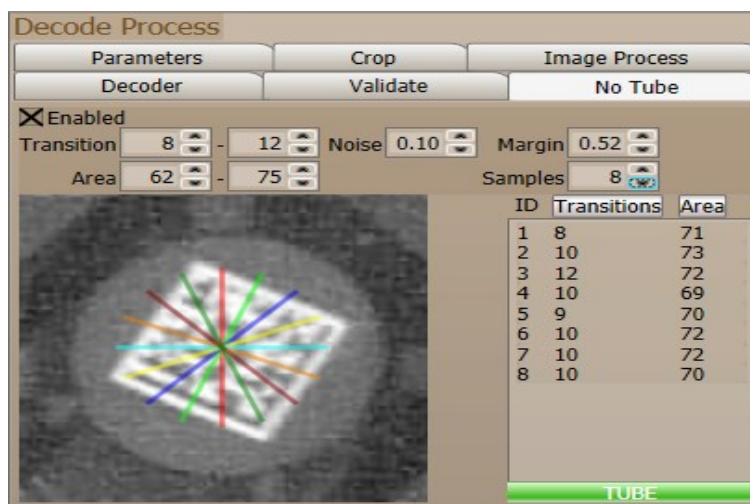

### Transitions

A transition occurs when the pixel value crosses the mean. A minimum and maximum value for the number of transitions on a sample line is required.

### Area

The modular area under the binaries for each sample line must fall between a minimum and maximum value to indicate that a matrix code may be present.

The reported Transitions and Area can be filtered numerically ascending or descending by selecting the Transition or Area white box.

The set value for each should allow the range to cover the min and max for each variable.


### Noise

Non-matrix code images often have many transitional lines scattered close to the mean (noise). These should be filtered out using the 'Noise Level' value, defined as a percentage of the range of pixel values in the sample.

### Margin

The width of a strip of pixels to be discarded around the edge of each cell image and is applied at decoding time. This should be set so the lines cross over the code and not the whole tube base.


## Decoder

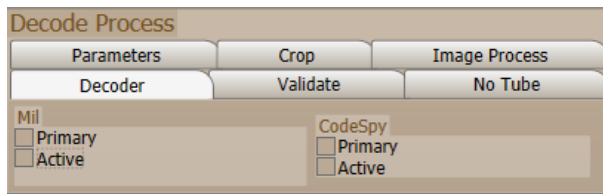

Option to select different decode methods

## Remote Connection via winsock

To communicate remotely to the device, add `-c` as a command line to the IntelliCode .exe in the IntelliCode folder.  
C:\Program Files (x86)\FluidX\IntelliCode  
Run the Application

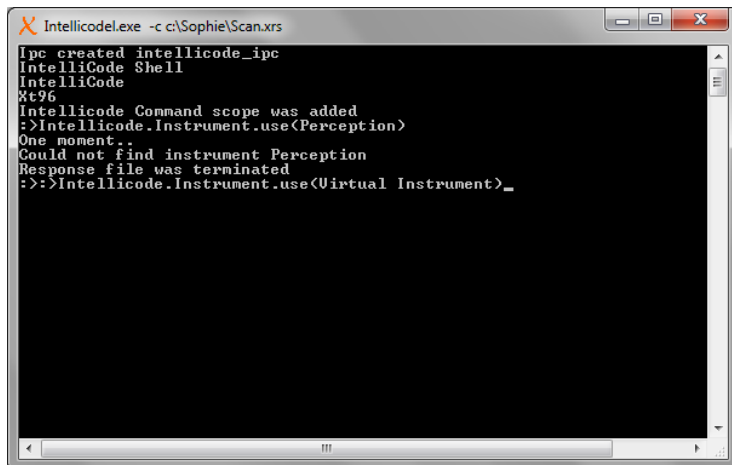

The software will try to open the last initiated device.

If not available the Instrument can be changed by using the following command

```
:>IntelliCode.Instrument.use(<instrument name>)
```

IntelliCode can auto fill text once the first few letters of a command are entered by pressing "Tab" on the keyboard.

For instance type "Int" Press "tab" and IntelliCode will automatically complete the rest of the text. "Intelli-

Code"

Selecting the "Up" key on the keyboard copies the last entry.

The commands are case and spelling sensitive. Please take care to enter information correctly or use the "Tab" shortcut.

**Info**= Retrieves information regarding Export name and types.

**list**= Retrieves list of available devices to run

**use**= Sets device

**Profile**=

