

CS

Submersible Fluorometer

User's Manual



September 19, 2011
P/N: 998-2300
Revision H


TURNER DESIGNS
845 W. Maude Ave.
Sunnyvale, CA 94085
Phone: (408) 749-0994
FAX: (408) 749-0998

C3 Submersible Fluorometer

Table of Contents

1. Introduction	
1.1 Introduction	4
1.2 Optical Sensors	4
1.3 Important Note Regarding the Use of the Fluorescein Sensor	5
2. Inspection and Setup	
2.1 Instrument Checklist	6
2.2 Optional Accessories	7
2.3 Identification	9
3. C-Soft Windows™ User Interface	
3.1 Software Installation and PC Requirements	10
3.1.1 Minimum PC Requirements	10
3.1.2 Software Installation	10
3.1.3 PC Software Interface	10
3.2 Software Operation	11
3.2.1 Settings Screen	11
3.2.2 Current Data and Calibration Screens	12
3.3 Calibration Procedures	
3.3.1 Direct Concentration Calibration	15
3.3.2 Raw Fluorescence Calibration – Blank Subtracted	16
3.3.3 Raw Fluorescence Calibration	17
3.3.4 Depth Calibration (Optional Accessory)	17
3.4 Log Setup Screen	18
3.5 Enabling Data Output Ready State	19
3.6 Downloading and File Management	20
4 Maintenance and Warranty	
4.1 Maintenance	21
4.1.1 Rinsing	21
4.1.2 Care for the Bulkhead Connector	21
4.2 Warranty Terms	21
4.3 Warranty Service	22
4.4 Out of Warranty Service	22
Appendices	
A C3 Submersible Fluorometer Specifications	24
B C3 Submersible Fluorometer Recommended Practices	25
C C3 Submersible Fluorometer Wiring and Cable Guide	26
D Linear Range, Quenching and Temperature Considerations	28
E C-FINS Installation Guide	29
F Installing the Battery Bracket	33
G Connecting the Submersible Battery Pack	34
H Installing the Mechanical Wiper	35
I Installing the Flow Cap	36
J Using the Secondary Solid Standard	38
K C-Series Analog Adapter	40
L How to set up Hyperterminal for Digital (ASCII) Data Capture	43

C3 Submersible Fluorometer

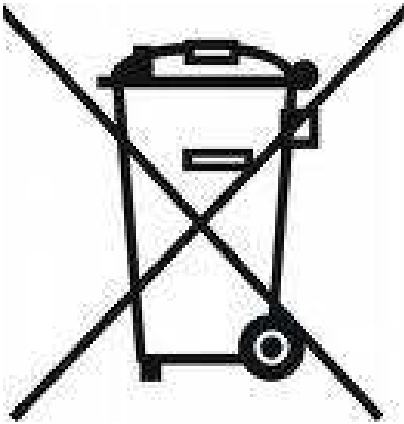

Equipment Specified as Electrical and Electronic Waste

C3 Submersible Fluorometer

1. Introduction

1.1 The C3 Submersible Fluorometer is manufactured according to users' optical specifications with one, two, or three optical sensors ranging in wavelengths from deep ultraviolet to infrared. Each optical sensor is designed with fixed excitation and emission filters. The C3 Submersible Fluorometer comes with a factory-installed temperature probe and can be configured with a pressure sensor for depth measurements up to 600 meters and/or a wiper to minimize biofouling. The C3 Submersible Fluorometer Windows™ based user interface allows for easy calibration of each sensor, digital data reporting, data logging and digital or analog export of data.

1.2 Optical configurations for Turner Designs' standard applications:

Application	Minimum Detection Limit	Dynamic Range	Excitation	Emission
<i>In Vivo</i> Chlorophyll <i>a</i>	0.025 µg/L	0-500 µg/L	430 nm	685/50 nm
Phycoerythrin (Marine Cyanobacteria)	150 cells/ml	0-150,000 cells/ml	540 nm	575/10 nm
Phycocyanin (Freshwater Cyanobacteria)	150 cells/ml	0-150,000 cells/ml	595 nm	> 645 nm
CDOM	0.4 ppb QS*	0-2500 ppb QS*	350 nm	430/30 nm
Turbidity	0.05 NTU	0-3000 NTU	850 nm	850 nm
Crude Oil	0.2 ppb PTSA**	0-2700 ppb PTSA**	350 nm	410-600 nm
Refined Fuels	2 ppb NS***	0-10,000 ppb NS***	260 nm	300-400 nm
Optical Brighteners	0.6 ppb PTSA**	0-15,000 ppb PTSA**	350 nm	445/15 nm
Fluorescein Dye (<i>See important note below</i>)	0.01 ppb	0-500 ppb	485 nm	540 nm
Rhodamine Dye	0.01 ppb	0-1000 ppb	550 nm	> 570 nm
PTSA Dye	0.01 ppb	0-650 ppb	365 nm	405/10 nm

* Quinine Sulfate in 0.05 Molar H₂SO₄

** 1,3,6,8-Pyrenetetrasulfonic Acid Tetrasodium Salt

*** 1,5-Naphthalene Disulfonic Disodium Salt

Custom optical configurations are available on request.

C3 Submersible Fluorometer

1.3 *Important Note Regarding Use of Fluorescein Optics:*

For C3's manufactured prior to 2/15/2011, the Fluorescein sensor should not be enabled for simultaneous data collection with Chlorophyll and/or Rhodamine sensors or it will produce unreliable results. For C3's manufactured after 2/15/2011 this interference issue was corrected. If you would like to send your C3 back to Turner Designs to resolve this interference issue, please contact sales@turnerdesigns.com

C3 Submersible Fluorometer

2. Inspection and Setup


2.1 Instrument Checklist

The C3 Submersible Fluorometer package (P/N: 2300-000) includes:

- ◆ C3 Submersible Fluorometer
- ◆ Factory-installed temperature sensor
- ◆ Interface cable for communicating with your PC and supplying power to the instrument
- ◆ CD including C-Soft software, user's manual, and quick start guide



12V Power Supply

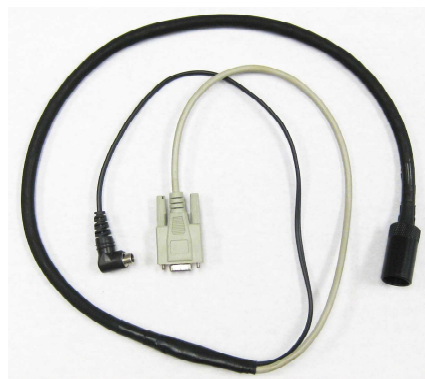

Computer Interface Cable

C3 Submersible Fluorometer

2.2 Optional Accessories

- ◆ Pressure sensor (P/N: 2300-360) *Factory-installed (Note: Can only be installed during the manufacturing of the unit.)*
- ◆ C3 Mechanical Wiper (P/N: 2300-450) *Factory-installed (Note: Can only be installed during the manufacturing of the unit.)*
- ◆ Replacement Wiper Brushes (P/N: 2300-460)
- ◆ Replacement Wiper Spacers (P/N: 2300-462)
- ◆ Rechargeable Submersible Battery Pack including charger (P/N: 2200-601)
- ◆ Battery Bracket (P/N: 2300-603) *(Note: Required for use with Battery Pack and Analog Output Adaptor.)*
- ◆ Flow Cap (P/N: 2300-700)
- ◆ Shade Cap (P/N: 2300-500) or C-ray Shade Cap (P/N 2300-502)
- ◆ Shade Cap Weight (P/N: 2300-510)
- ◆ C-ray Towed Deployment Body (P/N 2300-750)
- ◆ Extender Cables:
 - ◆ 10 meters (P/N: 105-2595)
 - ◆ 25 meters (P/N: 105-2596)
 - ◆ 50 meters (P/N: 105-2597)
 - ◆ Custom lengths available upon request
- ◆ Booster Kit (P/N: 2200-900) *(Note: Required for cable lengths greater than 10 meters.)*
- ◆ 24" Pigtail Interface Cable with Locking Sleeve (2200-170)
- ◆ 24" Integration Cable (P/N: 2200-160) or 5m Integration Cable (P/N: 2200-165) *(Note: Used for continuous data output and communication with the C3, NOT intended for uploading firmware to the C3.)*
- ◆ Secondary solid standard (P/N: 2300-900) for multiple sensors *(Note: Not available for turbidity or ultraviolet sensors.)*
- ◆ C-Series Analog Output Adapter (P/N: 2300-480)

Submersible Battery



Mechanical Wiper

C3 Submersible Fluorometer

Computer Integration Cable



C-ray Towed Deployment Body

Flow Cap



Shade Cap




Air Purge Slots

Battery Bracket



Booster Kit

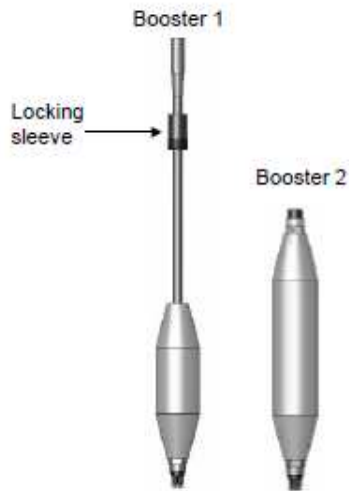

Solid Standard Insert



Secondary Solid Standard

C3 Submersible Fluorometer

2.3 Identification

Sensors are marked with numbers 1, 2, or 3 located adjacent to the optics on the face of the C3 Submersible Fluorometer. These numbers correspond to channels 1, 2, and 3 in C-Soft (C3 software). **Note: Channel names can be modified using C-Soft.**

Light emitting diode color table.

Application	Light Source Color
Chlorophyll a	Blue
Turbidity	Infra Red (No Color)
Phycocyanin	Yellow-Orange
Phycoerythrin	Green
Fluorescein	Blue <i>(See important note on page 5)</i>
Rhodamine	Green
CDOM	Ultra Violet (No Color)*
Refined Oils and Fuels	Deep Ultra Violet (No Color)*
Crude Oil	Ultra Violet (No Color)*
Optical Brighteners	Ultra Violet (No Color)*
PTSA	Ultra Violet (No Color)*

****Do not look directly at the optics. Ultraviolet light can be damaging to the eyes.***

C3 Submersible Fluorometer

3. C-Soft Windows™ User Interface

The C-Soft Windows™ based user interface allows for intuitive calibration, data logging, set up, and download functions.

3.1 Software Requirements and PC Installation

3.1.1 Minimum PC Requirements

- ◆ 133MHz microprocessor
- ◆ 600 X 800 VGA resolution
- ◆ Microsoft Windows™ 98
- ◆ 32 MB RAM

3.1.2 Software Installation

1. Exit all Windows™ programs
2. Insert the C-Soft CD (also located on the Turner Designs website on the Support page).
3. Open the C-Soft software folder and double click on the setup icon.
4. The software will automatically be installed in the PC's (C:/Program Files/C-Soft) folder unless user selects alternate location.
5. After restarting the PC double click on the "C-Soft" icon located on the desktop.



C-Soft Note: You may receive a warning message asking if you want to unblock messages, select "Unblock".


3.1.3 PC Software Interface

1. Connect the 9-pin serial end of the interface cable to the serial port on the back of the PC.

Note: If your computer only has USB ports we recommend purchasing a USB to serial adapter cable. <http://www.belkin.com/> search for Part Number: F5U409v1


2. Connect the 8-pin female end of the interface cable to the C3 Submersible Fluorometer's 8-pin male connector.
3. Connect the interface cable's 12V pin to the 12V power supply.
4. Plug the 12V power supply into a power source.
5. Wait for initialization process to complete.

C3 Submersible Fluorometer

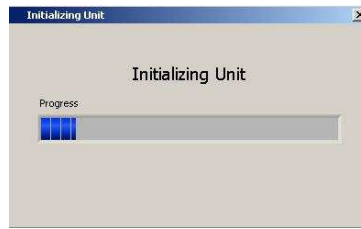

The connection icon, located on the upper right of the screen, will turn green if connected.

Note: If the connection icon does not turn green, disconnect and reconnect power.

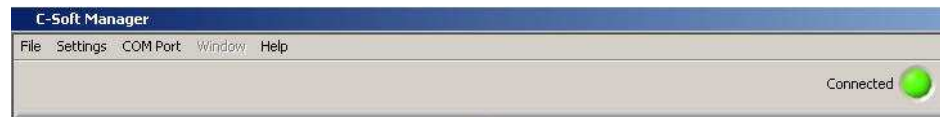


3.2 Software Operation

3.2.1 Settings Screen

The Settings Screen allows users to set the time and date, identify each optical sensor, update serial numbers, and enable/disable sensors.

- ◆ Users can manually set the date and time or choose to synchronize the C3 Submersible Fluorometer to the PC date and time.
- ◆ Each channel (1-3) in the Settings Screen has a location associated with it on the C3 Submersible Fluorometer that is numerically labeled (1-3), respectively. Channels 4-6 are designated for the C6 Multi-Sensor Platform and are disabled for the C3.
- ◆ Click on the box adjacent to the channel to enable channel operation. If there is no sensor installed users will be unable to select that channel.
- ◆ Channels can be renamed using the dropdown selections or by manually entering a name.
- ◆ The serial number boxes are designed for the C6 Multi-Sensor Platform. There is no serial number associated with the sensors in the C3.
- ◆ Settings will automatically be saved after exiting the Settings screen.

C3 Submersible Fluorometer

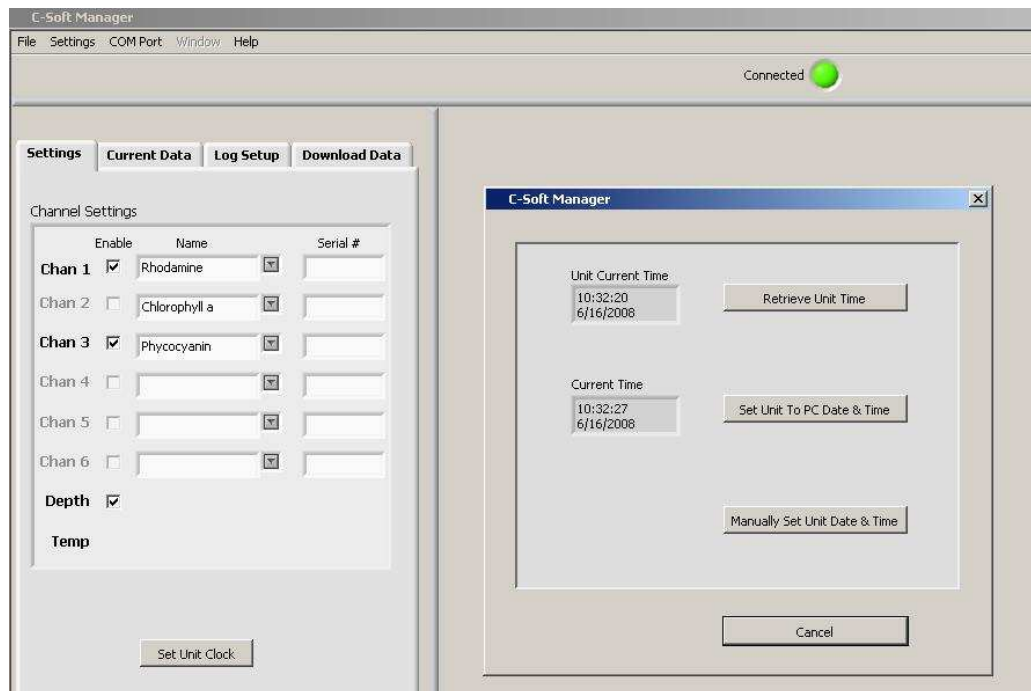

3.2.2 Current Data and Calibration Screen

The Current Data Screen allows users to view real-time data acquisition, in both tabular or waveform display, as well as conduct calibrations for each channel.


Current Data

- ◆ Channel names are displayed but cannot be changed in the Current Data screen.
- ◆ Real-time values are displayed next to each channel and in the adjacent Data Table.
- ◆ A calibration button is available for each channel and the depth sensor.
- ◆ Real-time data can be saved by selecting the “Start Data Capture” button. Users will be prompted to enter a file name and file location. Data files will be saved as .csv format.

Tabular Screen

- ◆ Real-time data including date, time, units and temperature are continuously scrolling on the tabular screen for enabled channels.
- ◆ Channel names will appear in the associated column headings.
- ◆ If units are not selected during calibration, a relative fluorescence units (RFU) label will appear.


C3 Submersible Fluorometer

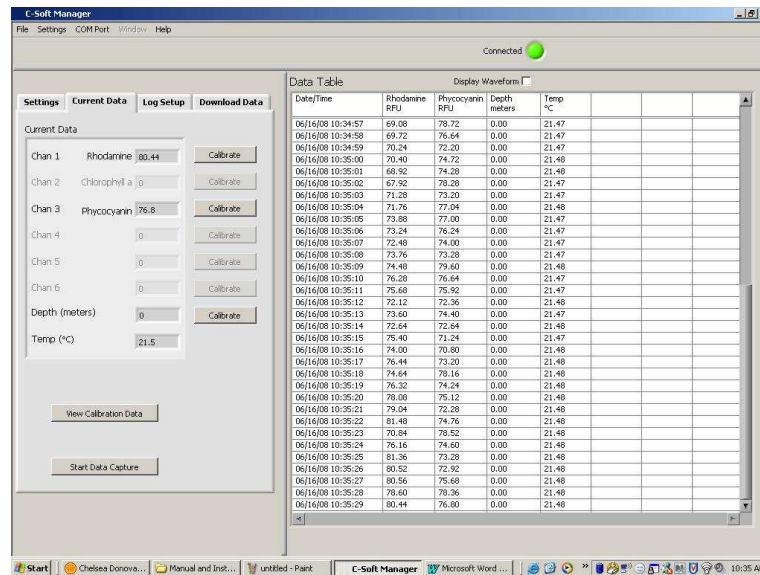

Waveform Screen

Users can graphically view real-time data for each sensor by clicking the “Display Waveform” box.

- ◆ Channel names and units will be displayed on the Y-axis.
- ◆ Time and date will be displayed on the X-axis.
- ◆ Clicking the “Autoscale” box will automatically scale graphs. Users may also rescale graphs by unchecking the “Autoscale” box and manually entering desired values for either Y or X axis ranges.
- ◆ Data can be saved during waveform mode by clicking “Start Data Capture”. Data files will be saved as .csv format.

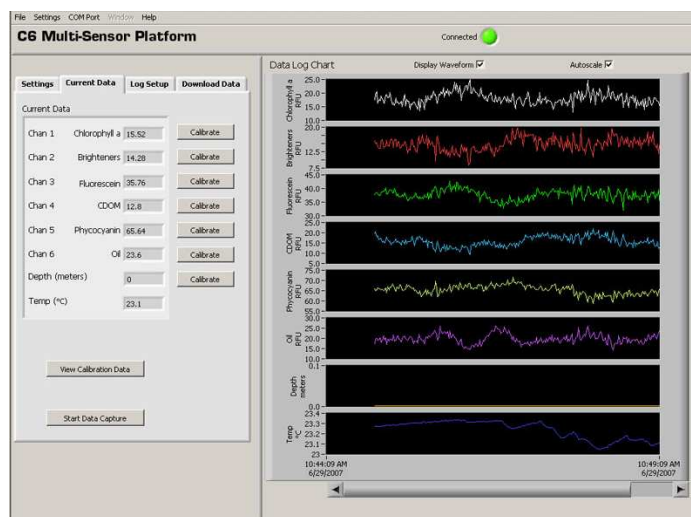

C3 Submersible Fluorometer

Calibration Screen

Users can calibrate sensors in either the direct concentration mode, raw fluorescence mode, or blank subtracted raw fluorescence mode.

Channel Calibration

Unit Label: Chan 1:Rhodamine Current RFU: 83.20

Step 1: Press "Set Blank" to set the blank value to the current RFU.

Blank RFU: x1 x10 x100

Step 2: Enter the value of your standard and select the units of measure.

Standard Value: Units:

Step 3: Press "Set Standard" to set the standard value to the current RFU.

Standard RFU:

Step 4: Optional: Click the temperature compensation box to activate.

Temp Compensation:

Step 5: After calibration is complete press "Save Calibration".

Direct Concentration Mode

Values reported in the Direct Concentration Mode are scaled to a predetermined standard value and blank subtracted. The equation used to calculate concentrations after calibrating to Direct Concentration Mode is:

$$\frac{\text{Standard Value (units)}}{\text{Standard RFU}} \times \text{Sample RFU} = \text{Concentration (units)}$$

Raw Fluorescence Blank Subtracted Mode

Values in the Raw Fluorescence Blank Subtracted Mode are noted as RFUB. It is important to note that these values are blank subtracted but are not scaled to a standard; they are **relative** values.

Raw Fluorescence Mode

Values in the Raw Fluorescence Mode are referred to as "Relative Fluorescence Units" (RFU). It is important to note that these values are not blank subtracted or scaled to a standard; they are **relative** values.

C3 Submersible Fluorometer

3.3 Calibration Procedure

3.3.1 Direct Concentration Calibration

Following steps 1-5 below will ensure values are reported as blank subtracted Concentrations in units of choice. See Appendix B for recommended practices.

Step 1: Blanking

- ◆ Place the C3 Submersible Fluorometer in a blank solution and ensure the optical head is free from air bubbles.
Note: A blank solution is a solution without the fluorophore of interest (i.e. de-ionized water, artificial water, or filtered water).
- ◆ Wait until “Current RFU” readings have stabilized then select the “Set Blank” button to set the blank value.
- ◆ Wait for the blank value to be measured and set for all 3 gains.

Step 2: Standard Value and Units

- ◆ Place the C3 Submersible Fluorometer in a standard solution or use a secondary solid standard (see Appendix D for more information on solid standards).
Note: UV and Turbidity sensors cannot be used with solid standards.
- ◆ Enter the known concentration value of the standard solution or solid standard in the “Standard Value” box. If the concentration is not known, enter an arbitrary value.
- ◆ Enter manually or select desired units from the dropdown menu.
Note: Do not choose or enter “RFUB” if Direct Concentration Mode is preferred.

Step 3: Set Standard

- ◆ When readings in “Current RFU” box have stabilized, select the “Set Standard” button.

Step 4: Optional Temperature Compensation

Temperature compensation is optional; it enables the C3 to compensate for changes in fluorescence due to varying temperatures as detected by the C3’s temperature sensor. Temperature compensation is available for Chlorophyll *a* and Rhodamine sensors only. Temperature is displayed in degree Celsius.

- ◆ Click on the “Temperature Compensation” box.
- ◆ Select either Chlorophyll *a* or Rhodamine from the dropdown menu.
- ◆ Data will now be corrected for temperature fluctuations.

C3 Submersible Fluorometer

The temperature compensation coefficients for each fluorophore are listed in the table below.

Fluorophore	Coefficient
Rhodamine	0.026 / °C Exponential
Chlorophyll <i>a</i>	0.014 / °C Linear

Step 5: Save Calibration

- ◆ After calibration is complete, select “Save Calibration”.
- ◆ Data are now displayed as concentration estimates.

3.3.2 Raw Fluorescence Calibration – Blank Subtracted

Following steps 1-5 below will ensure values are reported as Blank subtracted Relative Fluorescence Units (RFUB). See Appendix B for recommended practices.

Step 1: Blanking

- ◆ Place the C3 Submersible Fluorometer in a blank solution.

Note: A blank solution is a solution without the fluorophore of interest (i.e. de-ionized water, artificial water, or filtered water).

Note: Ensure the optical head is free of air bubbles.

- ◆ Wait until “Current RFU” readings have stabilized then select the “Set Blank” button to set the blank value.
- ◆ Wait for the blank value to be measured and set for all 3 gains.

Step 2: Selecting Units

- ◆ Select RFUB from the units’ dropdown menu.
Note: Do not enter a value in the standard value box.

Step 3: Skip and go to Step 4

Step 4: Optional Temperature Compensation – Follow same steps as the Direct Concentration Calibration procedure in Section 3.3.1.

Step 5: Save Calibration

- ◆ After calibration is complete select “Save Calibration”.
- ◆ Data are now displayed as blank corrected relative fluorescence units.

C3 Submersible Fluorometer


3.3.3 Raw Fluorescence Calibration

In the Current Data screen select the “calibrate” button next to a specific channel. At the bottom of the Calibration Screen select, “click to use Uncalibrated Raw Fluorescence Mode”. Data for that channel will be uncalibrated and reported as “Relative Fluorescence Units” (RFU).

3.3.4 Depth Calibration (Optional Accessory)

Depth is displayed in meters only. There are two calibration options for depth:

1. The first uses only the atmospheric pressure as an offset. This is suitable for any shallow application, including most fresh water applications.
2. The second also compensates for gravity variations with latitude, as designated from the UNESCO Technical Papers in Marine Science #44. This method assumes 0 °C and 35PSU salinity.

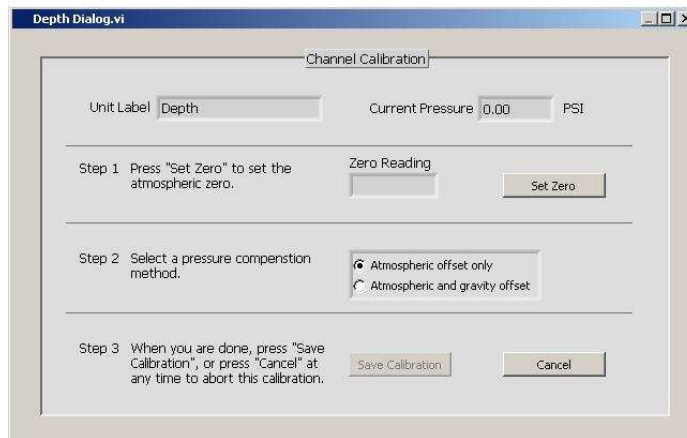

Atmospheric Offset

Step 1 (Method 1):

- ◆ Select the “Calibrate” button for depth in the “Current Data” screen.
- ◆ Pressure in PSI will be displayed in the “Current Pressure” window.
- ◆ After “Current Pressure” readings are stable at desired location select “Set Zero”.

Step 2 (Method 1):

- ◆ Select “Atmospheric offset only”.

Step 3 (Method 1):

- ◆ Select “Save Calibration” button.
- ◆ Depth values will now be scaled to the offset selected.

C3 Submersible Fluorometer

Atmospheric and Gravity Offset

Step 1 (Method 2):

- ◆ Select the “Calibrate” button for depth in the “Current Data” screen.
- ◆ Pressure in PSI will be displayed in the “Current Pressure” window.
- ◆ After “Current Pressure” readings are stable, at desired location, select “Set Zero”.

Step 2 (Method 2):

- ◆ Select “Atmospheric and gravity offset”.
- ◆ Enter the latitude.

Step 3 (Method 2):


- ◆ Select “Save Calibration” button.
- ◆ Depth values will now be scaled to the offset selected.

3.4 Log Setup Screen

The Log Setup screen allows the user to configure a logging session, set the mechanical wiper to its home position, and enable logging. The C3 Submersible Fluorometer can log just over 64,000 lines of data.

Enter a log start date (MM/DD/YYYY) and time (HH:MM:SS).

- ◆ Enter the sampling interval (HH:MM:SS). The minimum sampling interval is 1 second.

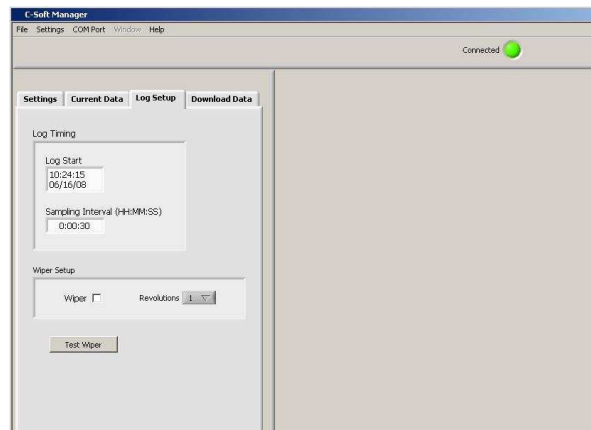


- ◆ If the mechanical wiper is installed click the “Wiper” box to enable.
- ◆ Select the number of revolutions from the dropdown box (1-10 revolutions). All sensors will be wiped at specified revolutions before each measurement.

Note: If users select revolutions that exceed sampling interval a warning message will be displayed.

Note: Do not click the Wiper Box to enable wiper if the wiper is not installed.

C3 Submersible Fluorometer

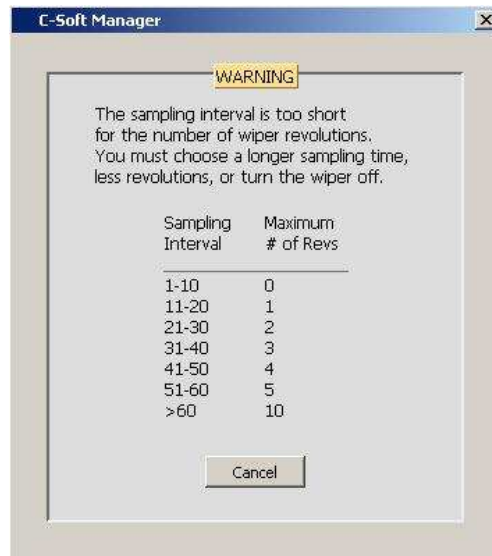

- ◆ After log set up is complete select “Enable Datalog” (see section 3.3 “Enabling Data Output Ready State”).
- ◆ A prompt will appear asking if user is ready to enable logging, select “Yes”.
Note: If data logging is set to start for some past time, there will be a minimum delay of 100 seconds before startup after power has been supplied. Otherwise wait until the data logging start time has been reached for data logging to begin.

WARNING! As soon as data logging is enabled users will be unable to access C-Soft settings. This should be the last step before the C3 Submersible Fluorometer is deployed.

3.5 Enabling Data Output Ready State

The Data Output Ready State must be enabled for the C3 Submersible Fluorometer in order to log data, capture real-time data output, and/or integrate the C3 with other instrumentation. There are five necessary steps to ensure the instrument is properly set up for data output:

1. Set the instrument’s clock to PC or local time in the Settings Tab.
2. Select the sensors you expect to use during sampling.
3. Calibrate selected sensors for either Direct Concentration Mode or Raw Fluorescence Units Blanked (RFUB) Mode (see section 3.3).
4. Set the start time, sampling, and wiping intervals in the Log Setup Tab.
Note: wiping interval is optional and depends on the sampling interval (see section 3.4).
5. Click the ‘Enable Datalog’ button followed by the ‘Yes’ button to enable the Data Output Ready State.

C3 Submersible Fluorometer

Note: Use the Battery Life Calculator available on the Turner Designs website C3 product page to calculate a conservative estimation of how many days you can expect to log data. Factors such as battery charge and water temperature will affect logging days.


A few seconds after the 'Yes' button has been clicked to enable the Data Output Ready State, the C-Soft program will automatically shut down and users must disconnect power and serial connections. The instrument is now in the Data Output Ready State and will remain in that state so long as C-Soft is not used to communicate with the instrument. **Note: As soon as C-Soft is used to communicate with the instrument, users will need to re-set the parameter settings in the Log Setup (section 3.4) and again enable datalog to put the instrument into the Data Output Ready State.**

When power is supplied to the instrument while it is in the Data Output Ready State, it will begin streaming data after a 100 second delay. If power is cycled (i.e. disconnected and reconnected) please allow 100 seconds for the data stream to begin.

3.6 Downloading and File Management

Users can download new and existing files saved in the C3 Submersible Fluorometer memory.

- ◆ Select Download Data.
Note: This step may take up to 10 minutes depending on file size.
- ◆ When download has completed choose the files to be saved by clicking the "Save" box.
- ◆ Users will be prompted to choose a folder location where data will be saved.
- ◆ Files will be saved in .csv format.
- ◆ Select "Clear Datalog" to erase existing data from memory.
- ◆ Users are able to graphically view downloaded .csv data files using C-Soft software if desired.

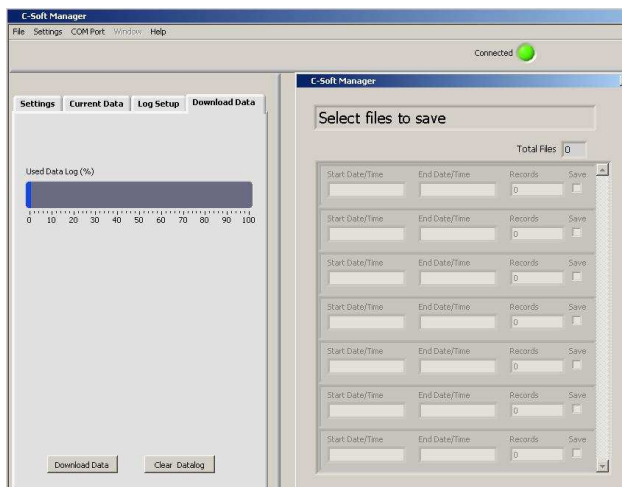

C3 Submersible Fluorometer

4. Maintenance and Warranty

4.1 Maintenance

4.1.1 Rinsing

The C3 Submersible Fluorometer should be rinsed or soaked in freshwater following each deployment.

4.1.2 Care for the bulkhead connector

Install the 8-pin female end plug to the bulkhead between uses. A light coat of Silicone spray should be used on the rubber of the male pins of the bulkhead to aid in sealing. The manufacturer recommends 3M™ Silicone Lubricant Spray or Loctite 8021 spray. **Note: You should avoid using silicone grease. Do NOT use WD-40, it will destroy the connectors.**

4.2 Warranty Terms

Turner Designs warrants the C3 and accessories to be free from defects in materials and workmanship under normal use and service for a period of 12 months from the date of shipment from Turner Designs with the following restrictions:

- Turner Designs is not responsible for replacing parts damaged by accident or neglect. Your instrument must be installed according to instructions in the User's Manual. Damage from corrosion is not covered. Damage caused by customer modification of the instrument is not covered.
- This warranty covers only Turner Designs products and is not extended to equipment used with our products. We are not responsible for accidental or consequential damages, except in those states where this limitation is not allowed. This warranty gives you specific legal rights and you may have other rights which vary from state to state.
- Damage incurred in shipping is not covered.

Extended warranties are available for purchase.
Contact sales@TurnerDesigns.com

C3 Submersible Fluorometer

4.3 Warranty Service

To obtain service during the warranty period, the owner shall take the following steps:

1. Write, email or call the Turner Designs Technical Support department and describe as precisely as possible the nature of the problem.

Phone: 1 (877) 316-8049

Email: support@turnerdesigns.com

2. Carry out any adjustments or tests as suggested by the Technical Support Department.
3. If proper performance is not obtained you will be issued a Return Materials Authorization number (RMA) to reference. Package the unit, write the RMA number on the outside of the shipping carton, and ship the instrument, prepaid, to Turner Designs. If the failure is covered under the warranty terms, the instrument will be repaired and returned free of charge, for all customers in the contiguous continental United States.

For customers outside of the contiguous continental United States who purchased equipment from one of our authorized distributors, contact the distributor. If you purchased directly, contact us. We will repair the instrument at no charge. Customer pays for shipping, duties and documentation to Turner Designs. Turner Designs pays for return shipment (custom duties, taxes and fees are the responsibility of the customer).


4.4 Out-of-Warranty Service

Follow steps for Warranty Service as listed above. If our Technical Support department can assist you by phone or correspondence, we will be glad to, at no charge. Repair service will be billed on a fixed price basis, plus any applicable duties and/or taxes. Shipment to Turner Designs should be prepaid. Your bill will include return shipment freight charges.

Address for Shipment:

Turner Designs, Inc.
845 W. Maude Ave.
Sunnyvale, CA 94085

C3 Submersible Fluorometer

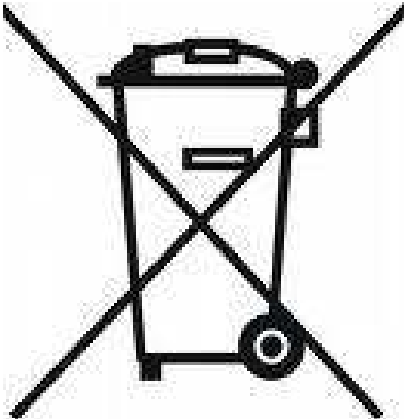

Equipment Specified as Electrical and Electronic Waste

Appendix A: C3 Submersible Fluorometer Specifications

C3 Submersible Fluorometer	
Weight in Air	1.64 kg (3.6 lbs)
Length	23 cm (9.1 in.)
Diameter	10 cm (3.9 in.)
Material	Delrin
Temperature	-2 to 50 °C
Temperature Accuracy	0.5 °C
Temperature Resolution	0.1
Depth	600 meters
External Power	8 to 30 VDC
Signal Output	Digital (ASCII) or Analog (<i>Note: analog requires accessory</i>)
Interface	RS232 (USB Adapter)
Maximum Sampling Rate	1 second
Nonvolatile Memory	64,000 Data Lines
Power Draw - Sleep Mode	< 1mAmp
Power Draw - Max	1 Watt

Submersible Battery Pack	
Battery	Lithium-Ion
Weight in Air	0.75 kg (1.65 lbs)
Weight in Water	0.22 kg (0.50 lbs)
Length	13.46 cm (5.3 in.)
Diameter	7 cm (2.75 in.)
Material	Delrin
Temperature	-2 to 50 °C
Depth	600 meters
Capacity	2,200 mAH
Voltage Output	14.8V nominal
Protection Features	Protects against: overcharge, over-discharge, overheating, short-circuiting

Appendix B: C3 Submersible Fluorometer Recommended Practices

Identification


Sensors are labeled 1, 2 or 3 adjacent to the optics on the face of the instrument. Each channel is factory set in the C3 Windows Software.

Note: The channel labels can be modified in the C-Soft software.

Recommended Measurement Practices

The following steps will improve the accuracy and repeatability of your measurements, especially at low concentration levels:

1. Use a dark plastic or glass container for your standards and/or samples. (**Note: Some plastics may fluoresce and might interfere with measurements.**)
2. If using a glass beaker, place it on a non-reflective surface, preferably black.
3. Ensure that the sensor is **>3 inches above the bottom** of the container.
4. Ensure that the sensor is in the center of the container, and has **>2 inches clearance** between the sensor and the inside surface of the container.
5. If using a glass beaker, check if the optical surface of the sensor is free of air bubbles. Remove air bubbles if any are observed. If using a dark plastic container, run the wiper or your finger over the optical sensors to ensure the sensor heads are free of air bubbles.

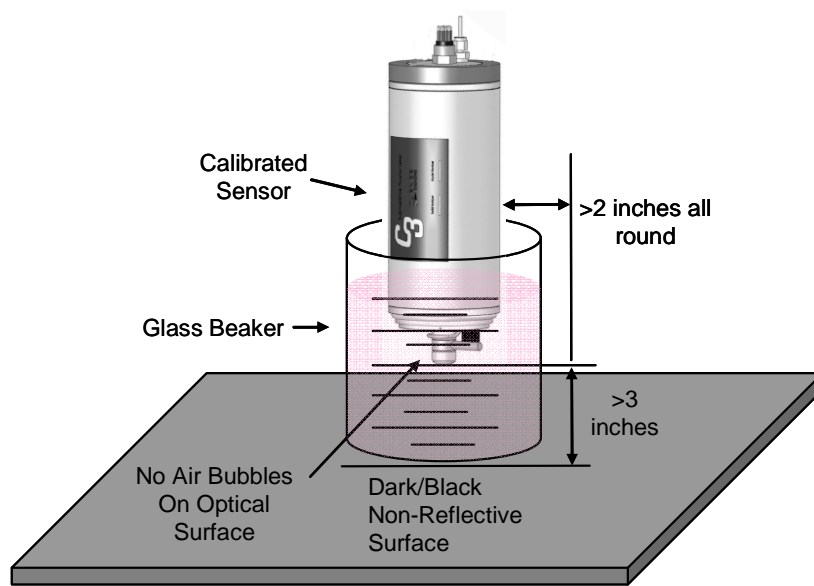


Appendix C: C3 Submersible Fluorometer Wiring and Cable Guide

The C3 Submersible Fluorometer outputs digital data in three formats: 1) Digital data that can only be read by C3 Submersible Fluorometer's software, 2) Digital data in ASCII format that can be read by larger multi parameter systems such as CTD's, and 3) Digital data that can be read and converted to analog data by the C-Series Analog Adapter.

An 8-pin Impulse cable provides two end connections: 1) a 9-pin RS232 serial port for connection to a PC or laptop computer and 2) a 12V port for supplying power to the unit.


C3 Submersible Fluorometer bulkhead and serial port connectors

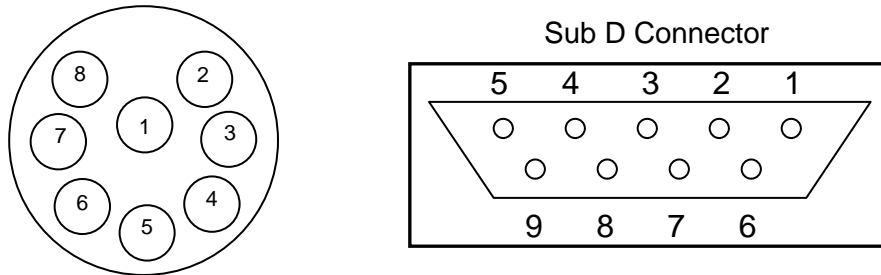

Wire Guide

Pin Out	Color	C3 Function	Corresponding Sub D Connector Pin
1	Black	V BATT (+)	Power Connector Cable – Center Pin (+)
2	White	V BATT (-)	Power Connector Cable – Housing (-)
3	Red	GND	PIN 5
4	Green	RX	PIN 2
5	Blue	TX	PIN 3
6	Brown	DTR & DSR	PIN 4, 6
7	Yellow	RTS & CTS	PIN 7, 8
8	Orange	N/A	N/A

**Power ground and V Batt (-) are not common.*

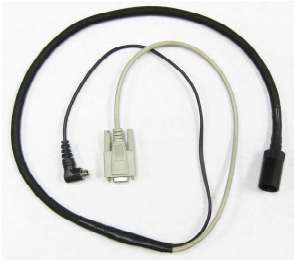

Computer Interface Cable

C3 Interface Cable (P/N 2200-150)

This cable is included as a standard accessory with the C3. It is used to program and calibrate the C3 and provide real-time data view. The cable length is approximately 2.5'.



Computer Integration Cable




yellow heat shrink

C3 Integration Cable (P/N 2200-160 and 2200-165) This cable is identified by the yellow band of heat shrink. The integration cable is used to connect the C3 to a PC or data logger for continuous data capture. It prevents interruptions in the C3 data transmission to the PC.



The cabled configuration allows for real-time data transmission. The C3 measurements are sent from the RS-232 cable to a PC where the fluorescence data can be viewed and recorded on a computer or a data logger that can record digital ASCII data. The data can also be saved on MS Excel from your C-Soft Software.

In order to provide greater length, extender cables P/N's 105-2595 (10 M), 105-2596(25M) or 105-2597(50 M) and custom cables up to (100M) can be installed along with the boosters (P/N 2200-900) here between the C3 and the interface cable to allow the signal and power to travel greater distances.

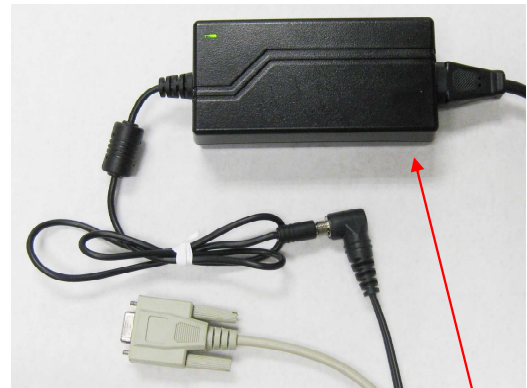

This is the +12VDC power supply P/N 159-0191. It is included as a standard accessory with the C3. The power supply connects to the +12 VDC connector on the interface or integration cable which will supply power to the C3.



This is where + DC power is connected.


This is the RS-232 connection to the PC

Appendix D: Linear Range, Quenching and Temperature Considerations

The linear range is the concentration range in which the C3 sensor output is directly proportional to the concentration of the signal. The linear range begins with the smallest detectable concentration and spans to an upper limit (concentration) that is dependent upon the properties of the material, filters used, and path length.

A non-linear relationship is seen at very high concentrations where the signal does not increase at a constant rate in comparison to the change in concentration (see figure below). At even higher concentrations, the signal will decrease even though the sample concentrations are continuing to increase. This effect is known as “signal quenching”.

Linearity can be checked by diluting a sample 1:1 or some other convenient ratio. If the sample is still in the linear range, the reading will decrease in direct proportion to the dilution. If the reading does not decrease in direct proportion to the dilution or if the reading increases, the sample is beyond the linear range.

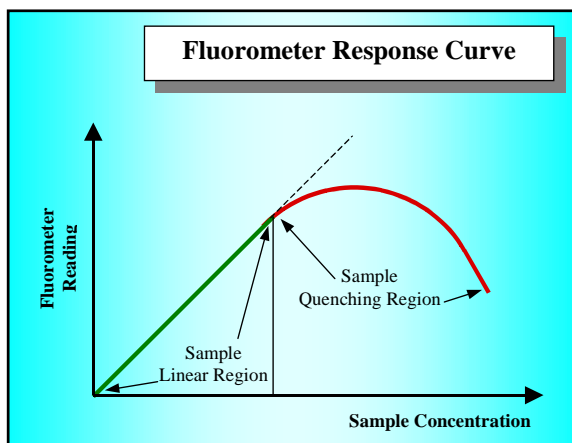

Graph showing Linear and Quenching Regions of the sample's response

Temperature Considerations

Fluorescence is temperature sensitive. As the temperature of the sample increases, the fluorescence decreases. For greatest accuracy, record the sample temperature and correct the sensor output for changes in temperature.

For further information on how temperature, light, water quality and the physiological state of the algal cells can all affect the measurement of chlorophyll *a*, please refer to the application section of Turner Designs' website.

Appendix E: CFINS Installation Guide

Requirements:

Turner Designs C3 Submersible Fluorometer with extender cable (boosters required for cable lengths >10M)
ArcGIS® Version 10
PC with two com (RS232 serial) ports or PCMCIA slot
GPS with com port (RS232) set to output NMEA and GPS serial cable
ArcMap C3 Extension


Steps:

Prepare C3 Fluorometer


- Run C-Soft and set up C3 for log sampling interval of 1 second (see Log Setup Section 3.4 of C3 User's Manual)
- Initiate datalog session as per instructions in Section 3.4 of manual. (C-Soft will automatically shut down)
- Power down and power up C3, it will take 90 seconds for data to start streaming

Prepare ArcMap


- If not already installed, install C3 extension in ArcMap (double click TD_C3.esriAddIn)
- Select both GPS Toolbar and Turner Designs C3 toolbar using *Customize/Toolbars*

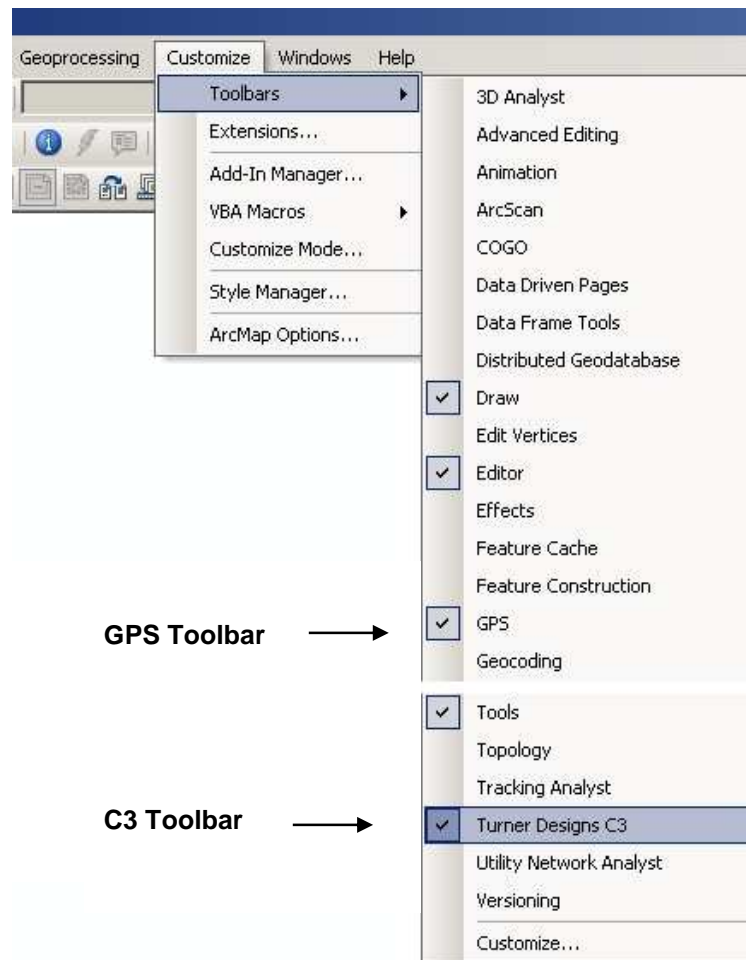


- Connect to GPS on GPS toolbar (See ArcMap help for information on this)

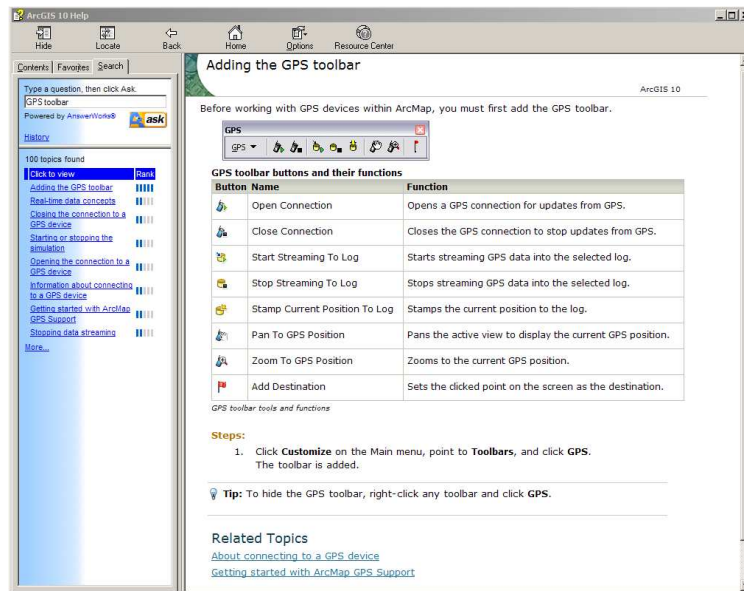

- Set up C3 Connection using C3 Toolbar (Setup C3)

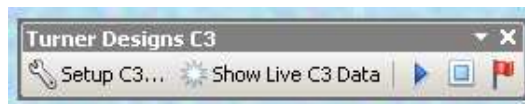

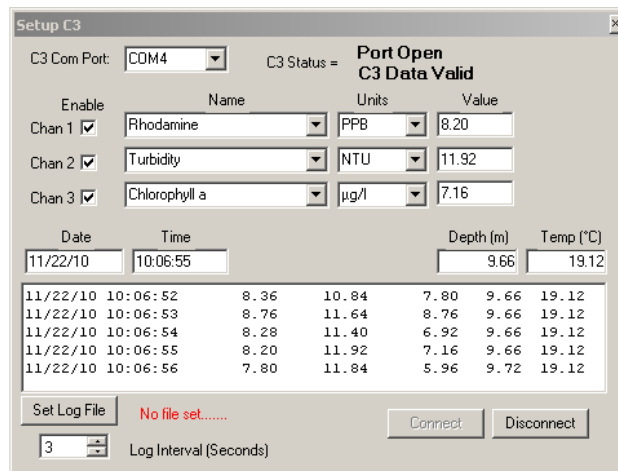


- Choose Com Port and Channel Settings then hit connect button. Data should start streaming in as shown below:



- Choose Shapefile to log to using Set Log File Button. (form can now be closed)

- Open Show Live C3 Data form which shows larger version of C3 data along with GPS data and Port/Logging Status.

- Start/stop logging to file using toolbar buttons

Button	Name	Function
	Log Data To File	Starts Logging to file based on Log interval
	Stop/Pause Logging to File	Stops/Pauses logging
	Flag Data Point ¹	Adds Flag to current data point and gets user notes on flag (e.g. "Bad reading C3 out of water" or "Start Transect 3")

¹ Note that flag only commits point once the OK button is pressed on user input. Long delays in writing notes may influence location of flagged point.

- Logged data has the following fields:

FieldName	Example	Notes
C3_Date	11/22/10	Current C3 Date
C3_Time	10:21:14	Current C3 Time
C3_DepthM	9.66	C3 Depth in Meters
C3_DegC	19.38	C3 Temperature in degrees Celsius
Chan1_Val	6.56	Channel 1 value
Chan1_Unit	PPB ²	Channel 1 Units (from C3 Setup Form. Not C-Soft)
Chan1_Type	Rhodamine	Channel 1 Name (from C3 Setup Form. Not C-Soft)
Chan2_Val	11.12	Channel 2 value
Chan2_Unit	NTU	Channel 2 Units (from C3 Setup Form. Not C-Soft)
Chan2_Type	Turbidity	Channel 2 Name (from C3 Setup Form. Not C-Soft)
Chan3_Val	7.68	Channel 3 value
Chan3_Unit	µg/l	Channel 3 Units (from C3 Setup Form. Not C-Soft)
Chan3_Type	Chlorophyll a	Channel 3 Name (from C3 Setup Form. Not C-Soft)
Lat_DD	-41.22914167	Latitude in Decimal degrees (Southern Latitudes negative)
Long_DD	173.0817033	Longitude in Degrees (Western Longitudes negative)
COG_True	17.58	Course over ground in degrees true
SOG	5.5556	Speed over ground in GPS Display Options Speed Units
SOG_Unit	Meters/Sec	GPS Speed over Ground Units
UTC_Date	21/11/2010	GPS date (in UTC)
UTC_Time	21:19:16	GPS Time (in UTC)
Flag	-1	Flag value (0 = no flag ; -1 = Flagged)
Flag_Notes	Start Transect 1	Flag Notes

² Channel names and units are logged as per the C3 setup form and not C-soft. These can be changed during a logging session but it is not recommended.


Appendix F: Installing the Battery Bracket

Introduction


The Battery Bracket (P/N: 2300-603) for the C3 Submersible Fluorometer was designed to provide support, allowing a direct connection between the battery and the C3 during self-contained deployment, thus eliminating the need for a cable connection. The Battery Bracket also provides balance and stability for the C3 Submersible Fluorometer and battery.

Installation

- 1) Use the extra plastic green support line provided with the Battery Bracket to remove the installed plastic green support line from the base cap of the C3 Submersible Fluorometer.

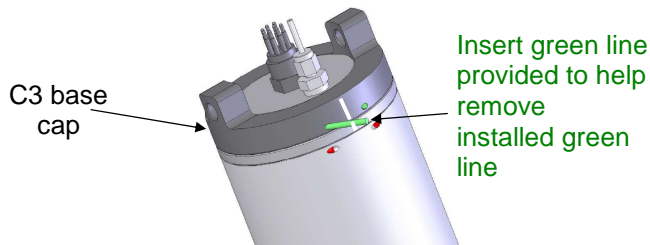


- 2) Pull the base cap off of the C3 Submersible Fluorometer.
- 3) Position the Battery Bracket with the C3 so the alignment strips line up with one another and the bracket's alignment pin fits into the C3's alignment hole.

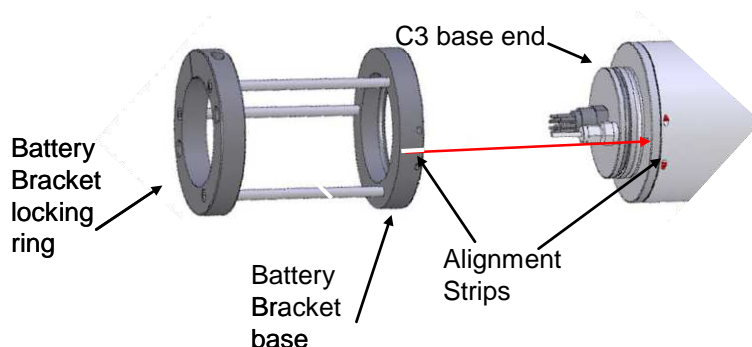

- 4) Secure the connection by feeding the Green Line through the holes on the Battery Bracket's base.



- 5) Slide the submersible battery pack through the bracket towards the male connector on the C3.
- 6) Mate the 8-pin male connector on the C3 with the female connector on the battery pack.
- 7) Secure the battery connection using the Allen Hex Key provided to tighten the Hex Screw located on the bracket's locking ring around the battery.

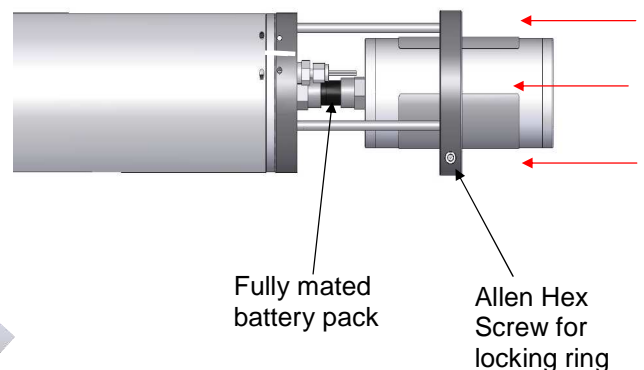

Appendix G: Connecting the Submersible Battery Pack

Introduction

The Submersible Battery Pack (P/N 2200-601) contains a rechargeable lithium-ion battery and battery pack charger. The battery pack is used together with the Internal Datalogger to provide self-contained use for the C6 Multi-Sensor Platform or C3 Submersible Fluorometer. The battery pack has a 600m depth rating and facilitates vertical profiling and moored applications.

Charging the Battery

- 1) Carefully align the battery's connector to the pins of the charger and push the battery straight onto the charger.
- 2) Plug the charger into a standard wall outlet and allow the battery to charge for 6 hours for a complete battery charge.
- 3) When finished, unplug the charger and remove the battery by pulling the battery straight out from the charger's connector.



General Precautions

Under no circumstances should you attempt to open the battery housing. Please contact Turner Designs if you have any problems or questions related to the battery pack.

Battery Life

The battery is designed, when fully charged, to exceed the Internal Data Logging capacity at the maximum sampling rate. For longer deployments at shorter sampling intervals, we recommend recharging the battery after the Internal Data Logger is full or every 3-4 months, whichever comes first.

Note: Use the Battery Life Calculator available on the Turner Designs website C3 product page to calculate a conservative estimation of how many days you can expect to log data. Factors such as battery charge and water temperature will affect logging days.

Specifications

Dimensions	5.3" x 2.75" (13.46cm x 7.0cm)
Weight in Air	1.65lbs. (0.75kg)
Weight in Water	0.50lbs (0.22kg)
Material	Delrin
Rated Depth	600m (1969 feet)
Voltage Output	14.8V nominal
Battery Type	Lithium-ion
Number of Cells	4
Capacity	2,200mAH
Charge Time*	6 hours
Connector	8 pin connector
Protection Features	Protects against: Overcharge, Over-Discharge, Overheating, Short-Circuiting

*Charge with Turner Designs charger only

Appendix H: Installing the Mechanical Wiper

Introduction


The Mechanical Wiper (P/N: 2300-450) for the C3 Submersible Fluorometer is designed to reduce bio-fouling during extended deployments.

Benefits

- Helps maintain consistent readings by reducing bio-fouling
- Replacement brushes available (P/N: 2300-460)


Installation

- 1) Connect the C3 to a computer that has C-Soft software and supply power.
- 2) Start the C-Soft software.
- 3) Click on the Log Data tab.
- 4) Click the Test Wiper button to bring the motor shaft to its home position.
- 5) Install O-ring into motor shaft groove. **Note: Do not lubricate O-ring**
- 6) Slide the wiper assembly onto the motor shaft with brushes facing towards the optical head and position the brushes between two sensors.

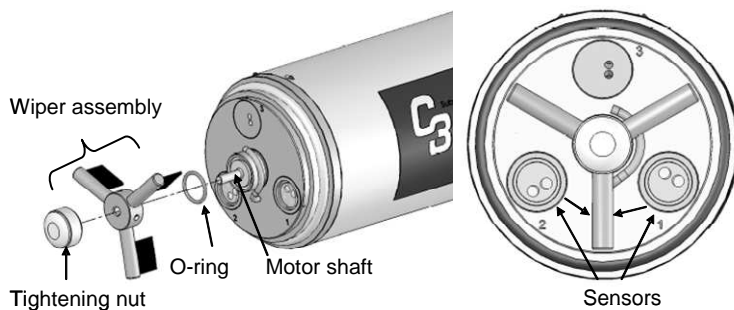


- 7) Hand-tighten nut to secure wiper assembly.

Note: do not use any tools to tighten the tightening nut.

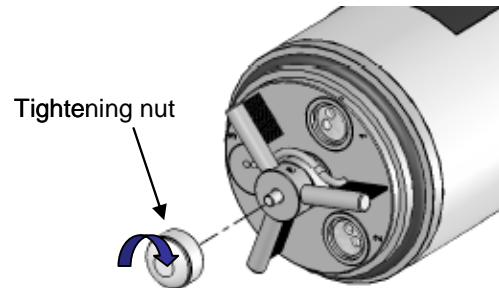

- 8) Click Test Wiper.
- 9) If the wiper does not come to rest in the desired position, unscrew the tightening nut, remove the wiper assembly, and repeat steps 6-8.

Replacement Brush Installation

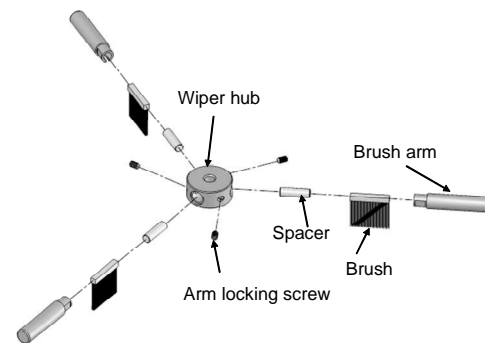

- 1) Using the Allen Wrench provided, loosen the arm locking screw.
- 2) Pull out the brush arm with brush.
- 3) Slide the brush towards the open end of the brush arm to remove the spacer and the brush.
- 4) Slide in a new brush. Slide in the spacer.
- 5) Position the brush arm onto the wiper assembly and tighten the arm locking screw.

Appendix I: Installing the Flow Cap

Introduction


The Flow Cap (P/N: 2300-700) for the C3 Submersible Fluorometer enables flow through mode. The Flow Cap will allow the C3 Submersible Fluorometer to be configured with other instruments used in flow through systems.

The C3 Flow Cap 1) provides elimination of other light sources that may interfere with sensor measurements, 2) protects sensors from damage during deployment or transport and 3) helps limit biofouling.

The Flow Cap can be installed on C3 Submersible Fluorometers with or without the mechanical wiper.

The Flow Cap can also be used as a calibration cup for C3 sensors using standard calibration solutions. See following page for calibration cup use.


Features

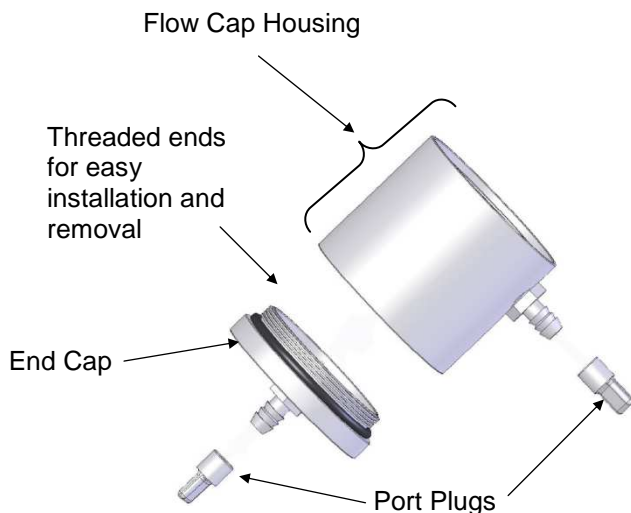

NOTE: The End Cap comes with an installed O-Ring that should not be removed or sealing will be compromised.


Installation Instructions

- 1) Turning clockwise, fully hand tighten the Flow Cap housing onto the C3 Submersible Fluorometer. The port located on the Flow Cap Housing should be closest to the optical head.

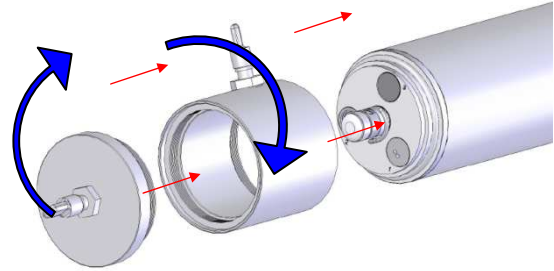

- 2) Turning clockwise, hand tighten the End Cap onto the Flow Cap Housing until it is fully tightened.
- 3) Flow Cap installation is complete.

NOTE: For optimal use, Turner Designs recommends positioning the C3 unit vertically when the Flow Cap is installed. This will purge any air from the system that might cause skewed readings.

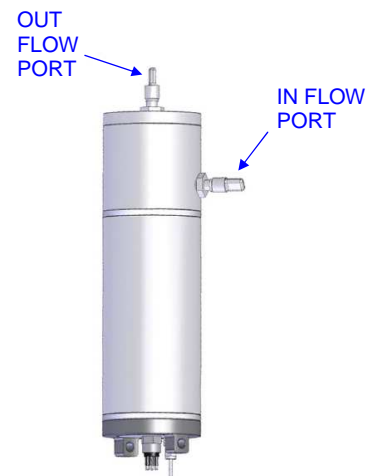


Calibration Cup Use

With the C3 Submersible Fluorometer's Flow Cap Housing installed:

- 1) Remove the End Cap.
- 2) Use a port plug to plug the port located on the Flow Cap Housing.
- 3) Position the unit vertically with optical head facing upward (see diagram to the right).

In this position the Flow Cap Housing can be used to hold standard solutions for calibrating the C3 sensors.

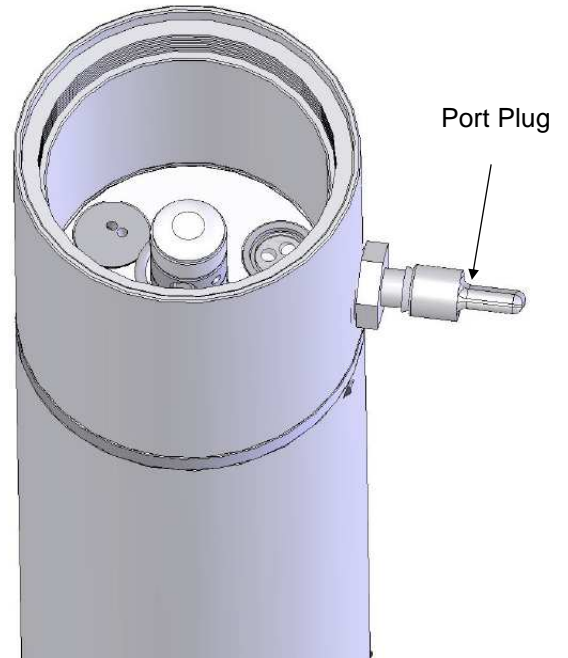

Specifications

Max. Pressure	100 psi
Length	10.1 cm (3.98 in.)
Diameter	10 cm (3.94 in.)
Weight	455 g (1.003 lbs.)
Max. Volume	400 ml
Material	Delrin

Appendix J: Using the Secondary Solid Standard

Introduction

The Solid Standard (P/N: 2300-900) for the C3 Submersible Fluorometer provides a very stable fluorescent signal which can be adjusted to:


- Match a specific sample
- Create a calibration correlation

Once a correlation is established between the Solid Standard and a liquid primary standard or typical sample it can be used for future calibrations. The Solid Standard can also be used as a stable reference point for checking signal stability and/or instrument drift.

Benefits


- Adjustable for a range of concentrations
- Can be used for quick and easy correlation.
- Can be used for multiple sensors
- Set screws ensure accurate positioning

Features

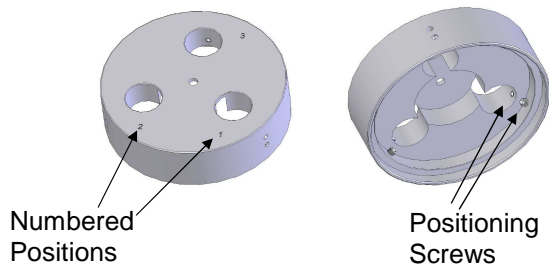


Installation and Use

- 1) To establish a correlation between a known concentration and its respective fluorescence output, immerse the sensor in a sample with a known concentration of your fluorophore.
- 2) Dry off the C3 Submersible Fluorometer's optics, attach the Solid Standard Cap.
- 3) Align Solid Standard Cap numbers with C3 sensor head numbers and snap on the Solid Standard Cap.

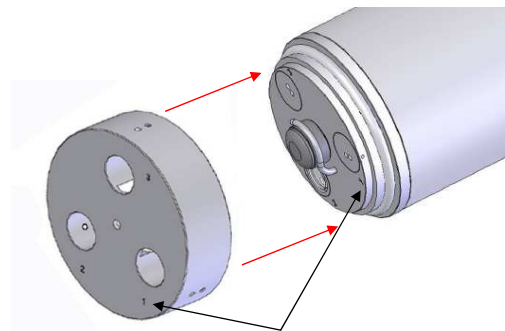

Solid Standard Cap numbers

- 4) Slightly rotate cap until it is set into position. **Note: You will feel/hear a click when cap is positioned.**
- 5) Insert up to 3 Solid Standards into the desired optical port position of the Solid Standard Cap.

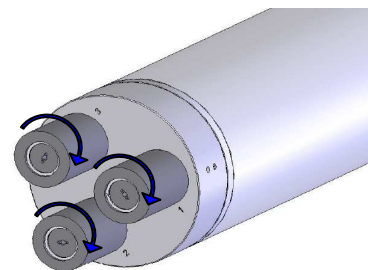


- 6) Rotate the Solid Standards until they set into position.

Note: You will feel/hear a click when the Solid Standard is positioned.

- 7) Adjust the Solid Standard to produce the same fluorescence output value from the sensor as in step 1.

To adjust each solid standard, use the green screwdriver to unscrew the locking nut as far as it will go.

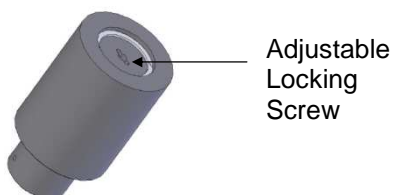

To change the signal level, insert the green screwdriver through the hole in the locking nut. Rotate it until it engages with the adjustment screw that is beneath the locking nut.

The signal can be adjusted to desired reading by turning the screw.

Once the desired reading has been obtained, the locking nut should be screwed down so that the adjustment screw is held firmly in place.

Note: The response of every solid standard is unique. Hence, a new correlation must be determined for every sensor.

- 8) Now your Solid Secondary Standard has been correlated to a known concentration and as long as it is no longer adjusted and is stored and used properly that

signal will remain a stable reference point.


- 9) For comprehensive information on fluorescence please refer to the application section of Turner Designs' website

Solid Standard Table

Application	Solid Standard
Chlorophyll a	Yes
Phycocyanin	Yes
Phycoerythrin	Yes
Rhodamine	Yes
Fluorescein	Yes
Solid standards are not available for Turbidity or Ultraviolet sensors	

Installation (C3 with wiper)

Remove the tightening nut and wiper arm from the C3.

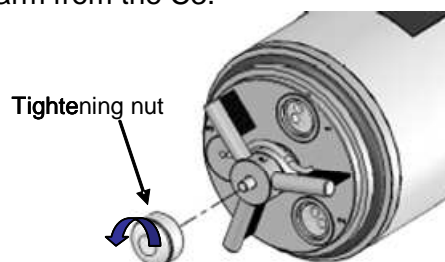

Now follow steps 1-4 of the installation instructions.

Ordering Instructions

Product	Part Number
Solid Standard Cap & 1 Solid Standard	2300-900
Individual Solid Standard	2300-901

Appendix K: C-Series Analog Adapter

Introduction

The C-Series Analog Adapter (P/N 2300-480) allows users to convert the standard digital output from the C3 Submersible Fluorometer to a 0 - 5 VDC analog signal. The Adapter can convert the signal from 1, 2, or 3 channels installed on a C3. The voltage output is proportional to the fluorescence intensity measured by the sensors. Modeled after the C3 Submersible Fluorometer's battery pack, the Analog Adapter is designed to connect directly to the C3. The C3 Battery Bracket (P/N 2300-603) is required to hold the Analog Adapter in place and secure it, preventing any break in communication or power. With the C-Series Analog Adapter users can integrate their C3 Submersible Fluorometers into any CTD or multi-parameter platform that will accept a 0 – 5 volt output signal.

Required Accessories

The following accessories are required for installation and operation of the Analog Adapter.

Battery Bracket (PN:2300-603) – Designed to hold the Analog Adapter in place and support it to avoid any break in communication or power ensuring continuous data output.

An Analog (Pigtail) Cable – We offer various lengths that are necessary for standard integrations.

- P/N: 2100-750 2 foot pigtail cable
- P/N: 2100-755 5 meter pigtail cable
- P/N: 2100-751 10 meter pigtail cable
- P/N: 2100-752 25 meter pigtail cable
- P/N: 2100-753 50 meter pigtail cable

C3 Submersible Fluorometer Setup

There are only a few steps required to set up the C3 Submersible Fluorometer for analog output. There are two methods that may be used for calibrating C3 channels/sensors. The first is a Raw Fluorescence Units Blanked Calibration (refer to section 3.3.2 in the C3's user manual) and the second is a Direct Concentration Calibration (refer to section 3.3.1 in the C3's user manual). Use the Raw Fluorescence Units Blanked (RFUB) calibration if you do not have a standard with known concentration or if you are unsure what numeric value to assign your standard.

NOTE: Repeat either one of the above calibrations for each channel that will be used for Analog Output.

Once calibration is complete, click the Log Setup tab to set the logging start time & logging interval. Click enable datalog, wait for C-Soft to automatically shutdown, and disconnect power. As soon as power is re-supplied to the C3 via the C-Series Analog Adapter it will automatically begin data output after 90 seconds, or pre-configured start time, at the set interval.

Analog Output

The C-Series Analog Adapter converts digital data to voltage to help users integrate our *in situ* fluorometers with almost any CTD, third party system, or data logger that accepts a 0-5 volt signal. Below are two ways to interpret voltage output based on the C3's setup.

Reading the Voltage Output for RFUB Calibrated C3's

The voltage output from the C3 Submersible Fluorometer via the C-Series Analog Adapter needs to be blank-corrected. Users can standardize the voltage signal using a standard of known concentration to convert RFUB to direct concentrations.

Reading the Voltage Output for Specific Range Calibrated C3's

The voltage output from the C3 Submersible Fluorometer via the C-Series Analog Adapter needs to be blank-corrected. The voltage output reflects direct concentrations re-scaled to 0-5 volts.

$$\text{Actual Concentration} = \frac{(\text{blank subtracted voltage})}{(5 \text{ volts})} \times (\text{concentration at 5 volts})$$

Example, if your measured voltage for a given sample is 2.567 volts and the blank voltage is 0.067 volts, your blank subtracted voltage is 2.5 volts.

$$[(2.5 \text{ volts}) / (5 \text{ volts})] = 0.5$$


If your range maximum equals 100 µg/L, then $(0.5 * 100 \text{ µg/L}) = 50 \text{ µg/L}$. Therefore, a sample that measures 2.567 volts = 50 µg/L.

Wiring Guide

The wiring guide below indicates the function of each pin for the Analog Adapter's 6-pin male end. To avoid damaging the Analog Adapter, please ensure the pins are only used for their intended purpose.

Pin	Wire Color	Function
1	Red	VBatt + (Positive)
2	Black	VBatt - (Negative)
3	White	Analog Out 1 (0-5 VDC)
4	Green	Analog Out 2 (0-5 VDC)
5	Blue	Analog Out 3 (0-5 VDC)
6	Brown	Analog GND


Analog Adapter's 6-pin male end

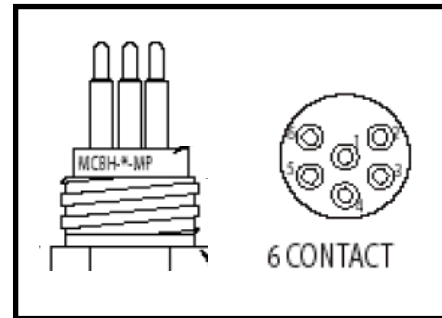


Appendix L: How to set up Hyperterminal for Digital (ASCII) Data Capture

1. Ensure C-Soft program is not running in the background. On an MS Windows computer, open the HyperTerminal program (Start→All Programs→Accessories→Communications→HyperTerminal)
2. The **Connection Description** screen will appear. Name the connection description. This file will save the communication parameters for the C3 Submersible Fluorometer and can be used in the future to establish communication with the instrument quickly.
3. The **Connect To** screen will appear. Choose the appropriate communications port in the **Connect Using** window. Click **OK**.
4. The **Port Settings** screen will appear. Choose the following port settings:

Bits per second: 9600
Data bits: 8
Parity: None
Stop bits: 1
Flow control: Hardware
Click **Apply** and **OK**

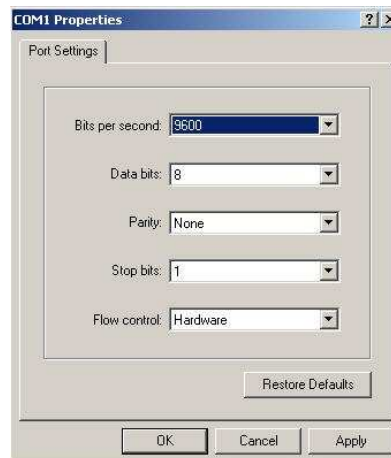

5. Next, the **HyperTerminal** window will appear. You are now ready to connect the C3 Submersible Fluorometer to a power source. Data will be displayed at set intervals in the HyperTerminal window after 100 seconds or at the specified log start time.

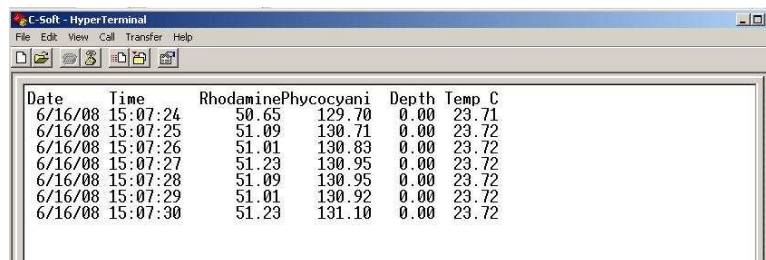

The image shows a screenshot of the 'C-Soft - HyperTerminal' window. The window displays a table of data with the following columns: Date, Time, Rhodamine, Phycocyanin, Depth, and Temp C. The data is as follows:

Date	Time	Rhodamine	Phycocyanin	Depth	Temp C
6/16/08	15:07:24	50.65	129.70	0.00	23.71
6/16/08	15:07:25	51.09	130.71	0.00	23.72
6/16/08	15:07:26	51.01	130.89	0.00	23.72
6/16/08	15:07:27	51.23	130.95	0.00	23.72
6/16/08	15:07:28	51.09	130.95	0.00	23.72
6/16/08	15:07:29	51.01	130.92	0.00	23.72
6/16/08	15:07:30	51.23	131.10	0.00	23.72

Most applications will integrate streaming data. Below are instructions on how to save streaming data on a PC.

To save data:

1. On the HyperTerminal window toolbar, choose Transfer and then Capture Text.



2. When you click on Capture Text, a window will appear asking you to **name** and **save** the **.txt** file to the location of interest.



3. Once you are finished, you will need to return to the HyperTerminal window toolbar, choose **Transfer** and then **Capture Text** and select **Stop**. This will stop data capture to the **.txt** file.

