# **ILS-LM Series**

# High-Performance Linear Motor Stages











USER'S MANUAL

# Warranty

Newport Corporation warrants this product to be free from defects in material and workmanship for a period of 1 year from the date of shipment. If found to be defective during the warranty period, the product will either be repaired or replaced at Newport's discretion.

To exercise this warranty, write or call your local Newport representative, or contact Newport headquarters in Irvine, California. You will be given prompt assistance and return instructions. Send the instrument, transportation prepaid, to the indicated service facility. Repairs will be made and the instrument returned, transportation prepaid. Repaired products are warranted for the balance of the original warranty period, or at least 90 days.

# **Limitation of Warranty**

This warranty does not apply to defects resulting from modification or misuse of any product or part.

# **CAUTION**

Warranty does not apply to damages resulting from:

- Incorrect usage:
  - Load on the stage greater than maximum specified load.
  - Carriage speed higher than specified speed.
  - Improper grounding.
    - ¬ Connectors must be properly secured.
    - ¬ When the load on the stage represents an electrical risk, it must be connected to ground.
  - Excessive or improper cantilever loads.
- Modification of the stage or any part thereof.

This warranty is in lieu of all other warranties, expressed or implied, including any implied warranty of merchantability or fitness for a particular use. Newport Corporation shall not be liable for any indirect, special, or consequential damages.

No part of this manual may be reproduced or copied without the prior written approval of Newport Corporation.

This manual has been provided for information only and product specifications are subject to change without notice. Any changes will be reflected in future printings.



**CAUTION** 

Please return equipment in the original (or equivalent) packing.


You will be responsible for damage incurred from inadequate packaging if the original packaging is not used.

# **Table of Contents**

Warra	ntyii
EC De	claration of Conformityv
Definit	tions and Symbolsvi
Warni	ngsvii
Cautio	onsix
1.0	— Introduction1
2.0	— <b>Description</b> 1
2.1	Design Details
3.0	— Characteristics
3.1	Definitions
3.2	Mechanical Specifications4
3.3	Hard Stoppers: Speed Limitation Versus Load4
3.4	Load Characteristics and Stiffness5
3.5	Example of Parameters "ScalingAcceleration" & "AccelerationLimit"5
3.6	Estimated Moving Mass for (M-)ILS-LM Stages
3.7	Stage Weights8
4.0	— Motor
4.1	Brushless Motor Characteristics8
4.2	Feedback Signal Position8
4.3	Pinouts9
	Motor Connector (SUB-D15M)9
	Encoder Connector (SUB-D15F)9
5.0	— Connection to Newport Controllers10
5.1	Warnings on Controllers10
5.2	Connection11
5.3	Cables11
6.0	— Connection to Non-Newport Electronics12
6.1	Connections
7.0	— Dimensions

9.0	— Maintenance	14
9.1	Maintenance	14
9.2	Repair	14
	Calibration	
Servic	e Form	15

# **EC Declaration of Conformity**

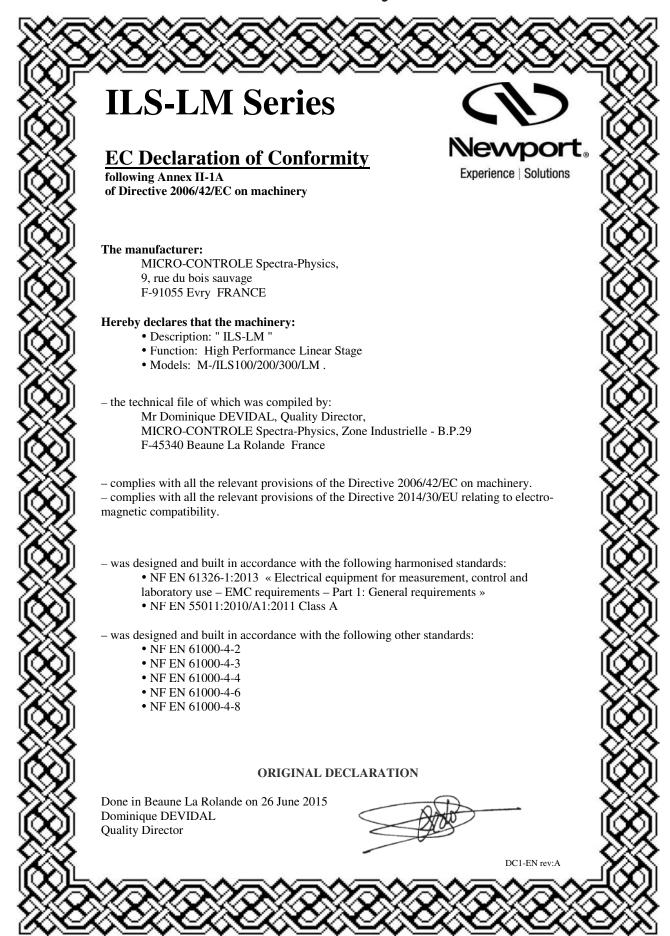

# **Definitions and Symbols**

The following terms and symbols are used in this documentation and also appear on the product where safety-related issues occur.

# **General Warning or Caution**



The exclamation symbol may appear in warning and caution tables in this document. This symbol designates an area where personal injury or damage to the equipment is possible.

The following are definitions of the Warnings, Cautions and Notes that may be used in this manual to call attention to important information regarding personal safety, safety and preservation of the equipment, or important tips.



#### **WARNING**

Warning indicates a potentially dangerous situation which can result in bodily harm or death.



## **CAUTION**

Caution indicates a potentially hazardous situation which can result in damage to product or equipment.

## **NOTE**

Note indicates additional information that must be considered by the user or operator.

# **European Union CE Mark**



The presence of the CE Mark on Newport Corporation equipment means that it has been designed, tested and certified as complying with all applicable European Union (CE) regulations and recommendations.

## **Warnings and Cautions**



# ATTENTION

This stage is a Class A device. In a residential environment, this device can cause electromagnetic interference. In this case, suitable measures must be taken by the user.

# Warnings



#### WARNING

The motion of objects of all types carries potential risks for operators. Ensure the protection of operators by prohibiting access to the dangerous area and by informing the personnel of the potential risks involved.



#### **WARNING**

The magnetic channel included in this device has the potential to disrupt pacemakers. Consequently, it is recommended that individuals maintain a distance of 1 meter or more from the stage as a precautionary measure.

#### WARNING

Do not use this stage when its motor is emitting smoke or is unusually hot to the touch or is emitting any unusual odor or noise or is in any other abnormal state.

Stop using the stage immediately, switch off the motor power and then disconnect the electronics power supply.

After checking that smoke is no longer being emitted contact your Newport service facility and request repairs. Never attempt to repair the stage yourself as this can be dangerous.

#### WARNING

Make sure that this stage is not exposed to moisture and that liquid does not get into the stage.



Nevertheless, if any liquid has entered the stage, switch off the motor power and then disconnect the electronics from power supply.

Contact your Newport service facility and request repairs.

#### **WARNING**

Do not insert or drop objects into this stage, this may cause an electric shock, or lock the drive.

Do not use this stage if any foreign objects have entered the stage. Switch off the motor power and then disconnect the electronics power supply.

Contact your Newport service facility for repairs.

# WARNING

Do not place this stage in unstable locations such as on a wobbly table or sloping surface, where it may fall or tip over and cause injury.

If this stage has been dropped or the case has been damaged, switch off the motor power and then disconnect the electronics power supply.

Contact your Newport service facility and request repairs.

# WARNING



Do not attempt to modify this stage; this may cause an electric shock or downgrade its performance.

# WARNING

Do not exceed the usable depth indicated on the mounting holes (see section "Dimensions"). Longer screws can damage the mechanics or cause a short-circuit.

# **WARNING**

Do not exceed speed and load limitations as specified in this manual.

# **Cautions**

#### **CAUTION**

Do not place this stage in a hostile environment such as X-Rays, hard UV,... or in any vacuum environment.

#### **CAUTION**

Do not place this stage in a location affected by dust, oil fumes, steam or high humidity. This may cause an electric shock.

#### **CAUTION**

Do not leave this stage in places subject to extremely high temperatures or low temperatures. This may cause an electric shock.

- Operating temperature: +10 to +35 °C
- Storage/Operating altitude: 1000 m
- Storage/Operating humidity: 85%
- Storage temperature: -10 to +40 °C (in its original packaging)



# **CAUTION**

Do not move this stage if its motor power is on.

Make sure that the cable to the electronics is disconnected before moving the stage. Failure to do so may damage the cable and cause an electrical shock.

# **CAUTION**

Be careful that the stage is not bumped when it is being carried. This may cause it to malfunction.

# **CAUTION**

When handling this stage, always unplug the equipment from the power source for safety.

# **CAUTION**

When the carriage is in its end-of-run position, it is strongly recommended not to go beyond this point as this may damage the stage mechanism.

# **CAUTION**

Contact your Newport service facility to request cleaning and specification control every year.

# High-Performance Linear Motor Stages ILS-LM Series

1.0

# Introduction

This manual provides operating instructions for the stage that you have purchased in the ILS-LM Series.



ILS300LM Stage.

## RECOMMENDATION

We recommend you read carefully the chapter "Connection to electronics" before using the ILS-LM stage.

2.0

# **Description**

The ILS-LM Series linear stages are designed for self-supporting applications with travel ranges from 100 mm to 300 mm. The stages feature a robust design with high performance but with low cost, making them cost-effective solutions for precision industrial applications such as semiconductor wafer inspection, micro-electronics test and assembly, pick and place, DNA sequencing, or laser machining,

The ILS-LM series linear motor stages are designed for self-supporting applications with travel ranges from 100 mm to 300 mm. The stages feature a robust design with high performance but with low cost, making them cost-effective solutions for precision industrial applications such as semiconductor wafer inspection, micro-electronics test and assembly, pick and place, DNA sequencing, or laser machining.

Based on industry proven technology of the ILS, the ILS-LM series utilizes an FEM-optimized extruded aluminum body that is extremely stiff, while minimizing the bending effect caused by the different thermal expansion coefficients of the aluminum body and the steel rails. The rails' position relative to the profile's neutral fiber minimize the effect due to bi-metal thermal expansion. The body's rigidity minimizes the deflection under load.

Unlike screw driven stages, the ILS-LM employs a centered, high efficiency 3-phase, synchronous ironless, linear motor as the driving element. This drive system is absolutely noise-free and has the advantage of higher speed, high acceleration and high system responsiveness without wear associated with motor brushes or drive screws. And because of the fully integrated linear motor, the ILS-LM is more than 150 mm shorter than a comparable screw driven stage. Thus, the ILS-LM is the optimum solution for space constrained applications that require high-throughput, high reliability, and ultra-quiet operation.

Recirculating ball bearing slides provide excellent payload capabilities and long life. The movement is smooth with low noise.

Precision position feedback is supplied by a very repeatable linear scale mounted in the stage. The encoder signals are interpolated by Newport's XPS motion controller with nanometer resolution for outstanding position sensitivity, repeatability, and stability. A home position is incorporated on the same scale, avoiding the use of additional electronics or mechanics for improved reliability and accuracy.

#### 2.1 Design Details

Base Material	Extruded Aluminum
Bearings	Recirculating bearings
Drive System	3-phase synchronous ironless linear motor (without Hall effect sensors)
Motor Initialization	Has to be done by the controller.
	The XPS controller uses an initialization routine to limit
	the stochastic motions of stages.
Motor Commutation	Done by the controller on encoder feedback
Feedback	Linear steel scale, 20 μm signal period, 1 V <sub>PP</sub>
Limit	Switches Optical
Home Switch	Optical, on encoder's fiducial track, located at center of travel
Controller Compatibility	ESP
Cable	5 m long cables included
MTBF	20,000 hours



# **NOTE**

This product complies with the RoHS directive (Restriction of Hazardous Substances).

#### **Characteristics**

#### 3.1 Definitions

Specifications of our products are established in reference to ISO 230 standard part II "Determination of accuracy and repeatability of positioning numerically controlled axes".

This standard gives the definition of position uncertainty which depends on the 3 following parameters:

## (Absolute) Accuracy

Difference between ideal position and real position.

# **On-Axis Accuracy**

Difference between ideal position and real position after the compensation of linear errors.

Linear errors include: cosine errors, inaccuracy of screw or linear scale pitch, angular deviation at the measuring point (Abbe error) and thermal expansion effects. All Newport motion electronics can compensate for linear errors.

The relation between absolute accuracy and on-axis accuracy is as follows:

Absolute Accuracy = On-Axis Accuracy + Correction Factor x Travel

### Repeatability

Ability of a system to achieve a commanded position over many attempts.

#### **Reversal Value (Hysteresis)**

Difference between actual position values obtained for a given target position when approached from opposite directions.

# **Minimum Incremental Motion (MIM or Sensitivity)**

The smallest increment of motion a device is capable of delivering consistently and reliably.

#### Resolution

The smallest increment that a motion device can theoretically move and/or detect. Resolution is not achievable, whereas MIM, is the real output of a motion system.

# Yaw, Pitch

Rotation of carriage around the Z axis (Yaw) or Y axis (Pitch), when it moves.

The testing of on-axis accuracy, repeatability, and reversal error are made systematically with test equipment in an air-conditioned room (20  $^{\pm 1}$  °C).

A linear cycle with 21 data points on the travel and 4 cycles in each direction gives a total of 164 points.

# **Guaranteed Specifications**

Guaranteed maximum performance values are verified per Newport's A167 metrology test procedure. For more information, please consult the metrology tutorial section in the Newport catalog or at **www.newport.com** 



# 3.2 Mechanical Specifications

Travel Range (mm)	100, 200, 300
Minimum Incremental Motion, Linear (μm)	0.01
Uni-directional Repeatability, Guaranteed (1) (μm)	0.45
Bi-directional Repeatability, Guaranteed (1) (µm)	0.7 or ±0.35
On-Axis Accuracy, Guaranteed (1) (2) (µm)	5 or ±2.5
Maximum Speed (mm/s) [See chapters 3.3 to 3.6]	500 (3)
Maximum Acceleration, No Load (m/s²) [See chapters 3.3 to 3.6]	20
Max. Force (Cont.) (N)	15
Max Force (Peak) 4 s (N)	30
Pitch, Yaw, guaranteed (1) (2) (4) (µrad)	300 or ±150
Load Capacity (N)	250



<sup>1)</sup> Shown are peak to peak, guaranteed specifications or ±half the value as sometimes shown. For the definition of typical specifications which are about 2X better than the guaranteed values, visit www.newport.com for the Motion Control Metrology Primer.

<sup>4)</sup> To obtain arcsec units, divide µrad value by 4.8.




# **CAUTION**

To reach specifications stated, stages must be fixed on a plane surface with a flatness of 5  $\mu m.$ 

# 3.3 Hard Stoppers: Speed Limitation Versus Load


ILS-LM stage uses electrical end-of-run and rubber bumper hard stoppers to stop as smoothly as possible the carriage at travel end. The overtravel allowed by the rubber bumpers is 3.5 mm.

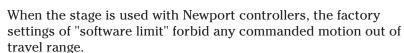

Nevertheless, for security issue, we recommend to follow the recommendations above to minimize risks of mechanical damage in case of failure or bad adjustment of parameters.

The maximum speed of the stage must be limited so that the rubber bumpers will always stop the carriage in 3.5 mm or less, to avoid any chock between carriage and stage body.

The graph opposite indicates the maximum speed that hard stopper can stop in 8 mm as a function of the load. This graph assumes that a correct wiring of the electrical end of runs will cut motor power before the contact with the rubber bumpers.

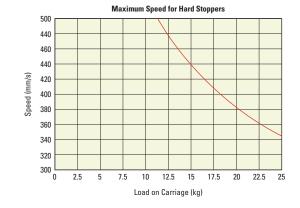


<sup>2)</sup> For a travel of 300 mm.

<sup>3)</sup> With a 44 V driver (such as the XPS-DRV02 board of the XPS controller).

#### 3.4 Load Characteristics and Stiffness

#### **Normal Load Capacity (Cz)**

Maximum load a stage can move while maintaining specifications.

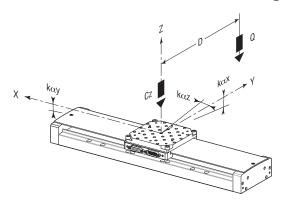

Cz, Normal centered load capacity	250 N
Kαx, Compliance in roll	15 μrad/Nm
Kαy, Compliance in pitch	10 μrad/Nm
Kαz, Compliance in yaw	10 μrad/Nm
Q, Off-center load (N)	0 ≤Cz ÷ (1 + D/60)
Where D = Cantilever distance (mm)	

# 3.5 Example of Parameters "ScalingAcceleration" & "AccelerationLimit"

This example is based on Newport XPS controller and XPS-DRV02 driver.

Two parameters have to be updated in the parameter file, according to the payload put onto the ILS-LM stage:

• The parameter "ScalingAcceleration" is used by the controller to scale the output voltage sent to the driver.

It indicates the theoretical maximum acceleration (friction not taken in account) of the ILS-LM carriage when the maximum voltage  $(10\ V)$  is applied to the driver.

• The parameter "AccelerationLimit" indicates the theoretical maximum acceleration that the controller will be able to require from the (M-)ILS-LM carriage. It correspond to the (M-)ILS-LM peak force.

For XPS controller and XPS-DRV02 driver, these parameters can be calculated using the following formulas, where the payload must be entered in kg:

ScalingAcceleration(Payload) = 
$$40305 \frac{\text{mm/s}^2}{\text{(Payload} + 0.9)}$$

**Examples:** ScalingAcceleration(0) =  $44800 \text{ mm/s}^2$ 

ScalingAcceleration(25) = 1556 mm/s<sup>2</sup>

$$AccelerationLimit(Payload) = \frac{ScalingAcceleration(Payload)}{1.1}$$

**Examples:** AccelerationLimit( $\mathbf{0}$ ) = 40.727 mm/s<sup>2</sup> AccelerationLimit( $\mathbf{25}$ ) = 1712 mm/s<sup>2</sup>

These formulas are explained below:

#### Inputs from the system

• Motor:

Motor force constant: MotorForceConstant = 11.4 N/Amp. rms

- Motor thermal resistance: MotorRth = 1.8  $\frac{K}{Watt}$ 

- Motor constant at 300 K: MotorK =  $18.3 \frac{\text{Newton}^2}{\text{Watt}}$ 

#### • Driver:

XPS-DRV02 driver is using sine/cosine commutation on its input. Its maximum current of 5 A is obtained for a 10 V input, so:

MaxDriverCurrent = 5 Amp.

TransImpedenceDriver = 0.5 Amp./V

#### Mechanics:

ILS-LM carriage mass: Mcar = 0.4 kgInterface plate mass: Mint = 0.5 kg

# • System:

ILS-LM rms force is defined to keep motor heating below 25 °C:

$$ILS\_RmsForce = \sqrt{\frac{25 \cdot MotorK}{MotorRth}} \\ ILS\_RmsForce = 15 \ Newtons$$

ILS-LM peak force is defined to allow two time the rms force. This value is lower than the theoretical peak force of the motor:

The rms current limitation of the driver is set according to (M-)ILS-LM rms force:

The peak current limitation of the driver is set according to ILS-LM peak force:

$$\label{eq:decomposition} \begin{aligned} & \text{DriverMaximumPeakCurrent} = min \bigg( \frac{ILS\_PeakForce}{MotorForceConstant} \sqrt{2} \cdot 1.05 \text{ , MaxDriverConstant} \bigg) \\ & \text{DriverMaximumPeakCurrent} = 3.9 \text{ A} \end{aligned}$$

The integration time of the driver  $I^2t$  limitation is set to allow to apply the peak current for 4 seconds. This is obtained with an integration time of 10 s, which is lower than the motor thermal time constant.

DriverRMSIntegrationTime = 10 s

# Inputs from the user

Load on the carriage: Load = 0 kg

#### **Calculation**

# MovingMass:

MovingMass = Mcar + Mint + Load

### • ScalingAcceleration:

$$ScalingAcceleration = \frac{\text{MotorForceConstant} \cdot 10 \text{ Volts} \cdot \text{TransImpedenceDriver}}{\sqrt{2} \cdot \text{MovingMass}}$$
 
$$ScalingAcceleration = \frac{\text{44800 mm/s}^2}{\text{Voltage Move MovingMass}}$$

# • LimitAcceleration:


LimitAcceleration = ScalingAcceleration • DriverMaximumPeakCurrent • 1.05
LimitAcceleration = 23670 mm/s<sup>2</sup>

# 3.6 Estimated Moving Mass for (M-)ILS-LM Stages





# **CAUTION**


ILS-LM stages with linear motor are sensitive to the load variation and its stiffness. The example below shows the behavior of the displacement in accordance with the underestimated/overestimated moving mass.

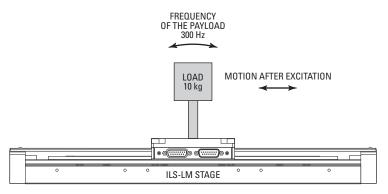

Driving in force requires a good adjustment of the controller/driver theoretical acceleration parameter (ScalingAcceleration) which depends on the moving mass.

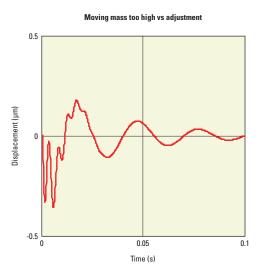

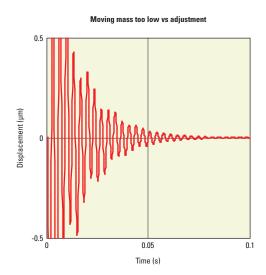

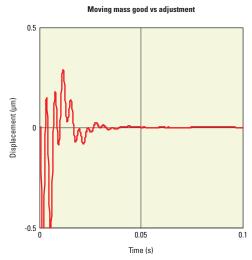

# 3.7 Stage Weights

Weights indicated into the below table are average values for stages with a typical drive unit installed.

The stage weight below does not include the cables.

	Weight [lb (kg)]
ILS100LM	6.4 (2.9)
ILS200LM	8.2 (3.7)
ILS300LM	9.9 (4.5)


4.0

# **Motor**

# 4.1 Brushless Motor Characteristics

Motor	Motor Constant	Magnet Pitch	Nominal Voltage	Max. RMS Current	Max. Peak Current	Resistance per Phase	Inductance per Phase
	$(N^2/W)$	(mm)	(V)	(A)	(A)	$(\Omega)$	(mH)
UC6	18.3	16.5	48	1.3	3.9	2.4	0.38


# 4.2 Feedback Signal Position

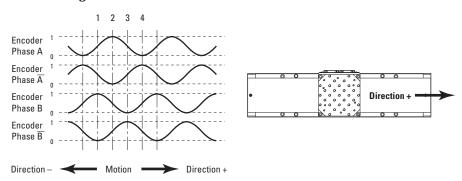

# 4.3 Pinouts


SUB-D connections for ILS-LM stages are given in the following tables:

# **Motor Connector (SUB-D15M)**

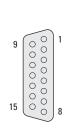

1	Shield Ground
2	V Motor
3	Mechanical Zero
4	- End-of-Run
5	Shield Ground
6	N.C.
7	N.C.
8	N.C.
9	W Motor
10	U Motor
11	+ End-of-Run
12	+5 V
13	N.C.
14	N.C.
15	N.C.

# **Encoder Connector (SUB-D15F)**

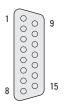

1	Encoder Phase A
2	Shield Ground
3	Encoder Phase B
4	Encoder Power: +5 V
5	N.C.
6	N.C.
7	Index Pulse /I
8	N.C.
9	Encoder Phase /A
10	N.C.
11	Encoder Phase /B
12	N.C.
13	N.C.
14	Index Pulse I
15	N.C.

# **Connection to Newport Controllers**

# 5.1 Warnings on Controllers

Controllers are intended for use by qualified personnel who recognize shock hazards and are familiar with safety precautions required to avoid possible injury. Read the controller user's manual carefully before operating the instrument and pay attention to all written warnings and cautions.

#### **WARNING**

Disconnect the power plug under the following circumstances:

- If the power cord or any attached cables are frayed or damaged in any way.
- If the power plug is damaged in any way.
- If the unit is exposed to rain, excessive moisture, or liquids are spilled on the unit.
- If the unit has been dropped or the case is damaged.
- If you suspect service or repair is required.
- Whenever you clean the electronics unit.

#### **CAUTION**

To protect the unit from damage, be sure to:

- Keep all air vents free of dirt and dust.
- Keep all liquids away from the unit.
- Do not expose the unit to excessive moisture (85% humidity).



# Read this manual before using the unit for the first time.

#### **WARNING**

All attachment plug receptacles in the vicinity of this unit are to be of the grounding type and properly polarized.

Contact your electrician to check your receptacles.

# WARNING

This product is equipped with a 3-wire grounding type plug.

Any interruption of the grounding connection can create an electric shock hazard.

If you are unable to insert the plug into your wall plug receptacle, contact your electrician to perform the necessary alterations to ensure that the green (green-yellow) wire is attached to earth ground.

# WARNING

This product operates with voltages that can be lethal.

Pushing objects of any kind into cabinet slots or holes, or spilling any liquid on the product, may touch hazardous voltage points or short out parts.

# 5.2 Connection

On each stage is represented a label which indicates its name and its serial number.



#### WARNING

Always turn the controller's power OFF before connecting to a stage.

Stages may be connected to the rear panel motor connectors any time prior to power-up with the supplied cable assemblies.

# **NOTE**



These stages are ESP compatible. Enhanced System Performance is Newport's exclusive technology that enables Newport ESP motion controllers to recognize the connected Newport ESP stage and upload the stage parameters. This ensures that the user can operate the motion system quickly and safely.

## 5.3 Cables

ILS-LM stages are delivered with a set of two 5-meter cables that can be directly connected to the XPS Newport controller.

# **WARNING**

ILS-LM Series translation stages can only operate with cable lengths of 5 m max.



## **WARNING**

These cables are shielded. For correct operation, make sure to lock connectors (ground continuity provided by cables).

## **WARNING**

Keep the cables at a safe distance from other electrical cables in your environment to avoid potential cross talk.

# 6.0

# **Connection to Non-Newport Electronics**

#### 6.1 Connections

# **WARNING**


Newport is not responsible for malfunction or damage to an ILS-LM stage when it is used with non-Newport controllers.


# **WARNING**

Newport guarantees "( "compliance of the ILS-LM stages only if they are used with Newport cables and controllers.

Nevertheless, the figure below shows the wiring when an ILS-LM stage is used with non-Newport controllers.




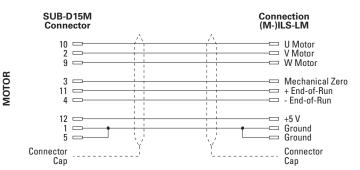

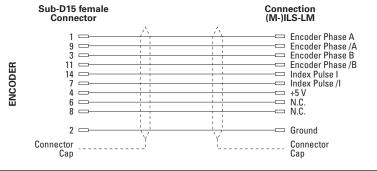

# 7.0

# **Dimensions**

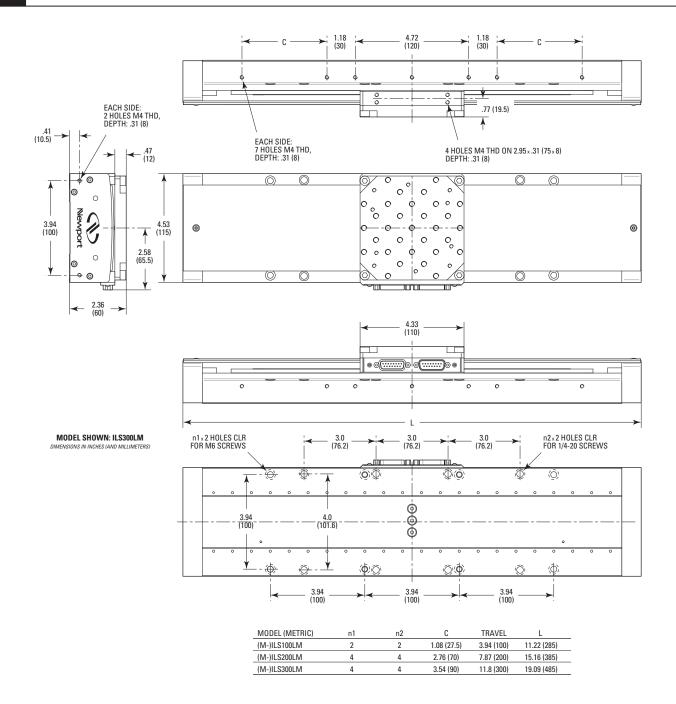

# 9.0

#### **Maintenance**

#### RECOMMENDATION

It is recommended to contact our After Sales Service which will know to define the appropriate maintenance for your application.

#### 9.1 Maintenance

The ILS-LM stage requires no particular maintenance. Nevertheless, this is a precision mechanical device that must be kept and operated with caution.

#### **PRECAUTIONS**

The ILS-LM stage must be used or stocked in a clean environment, without dust, humidity, solvents or other substances.

#### RECOMMENDATION

It is recommended to return your ILS-LM stage to Newport's After Sales Service after every 2000 hours of use for lubrication.

If your stage is mounted on a workstation and cannot be easily removed, please contact Newport's After Sales Service for further instructions.

# 9.2 Repair



## **CAUTION**

Never attempt to disassemble a component of the stage that has not been covered in this manual.

To disassemble a non specified component can cause a malfunction of the stage.

If you observe a malfunction in your stage, please contact us immediately to arrange for a repair.



# **CAUTION**

Any attempt to disassemble or repair a stage without prior authorization will void your warranty.

### 9.3 Calibration



## **CAUTION**

It is recommended to return your ILS-LM stage to Newport once a year for recalibration to its original specifications.

# **Service Form**

Name:	Return authorization #:	
Company:	(Please obtain prior to return of item)	
Address:		
Country:		
P.O. Number:		
Item(s) Being Returned:		
Model #:	Serial #:	
Description:		
	ny specific problems):	

**Your Local Representative** 

Tel.: \_\_\_\_\_

Fax: \_\_



# **Nevvport**®

# Experience | Solutions

# Visit Newport Online at: www.newport.com

# North America & Asia

Newport Corporation 1791 Deere Ave. Irvine, CA 92606, USA

# **Sales**

Tel.: (800) 222-6440

e-mail: sales@newport.com

# **Technical Support**

Tel.: (800) 222-6440

e-mail: tech@newport.com

# Service, RMAs & Returns

Tel.: (800) 222-6440

e-mail: service@newport.com

# **Europe**

MICRO-CONTROLE Spectra-Physics S.A.S

9, rue du Bois Sauvage 91055 Évry CEDEX France

# **Sales & Technical Support**

Tel.: +33 (0)1.60.91.68.68

e-mail: france@newport.com

# **Service & Returns**

Tel.: +33 (0)2.38.40.51.55