

User Manual

Sartorius Cubis Series

Electronic Semimicro, Micro, Precision and Analytical Balances MSA Models



Contents

Notes on Using this Manual	Weighing in Perc	ent	102
Safety Instructions	Timer-controlled	Functions	105
Intended Use	Totalizing		107
Overview of Models	DKD Measuremen	nt Uncertainty	109
Getting Started 8	Second Tare Men	nory (Preset Tare)	111
Transporting the Balance 10	Parts Counting		113
Installation Instructions 1	Checkweighing		117
Power Connection	Importing/Export	ting Data	121
Anti-theft Locking Device	Calibration and A	Adjustment	124
Below-Balance Weighing	Calibration/Adj	justment Using Internal Check Weight.	124
Using the Balance	Calibration/Adj	justment Using External Check Weight.	125
Turning the Balance On/Off	User Managemen	nt	127
Operating Design: Q-Guide	User Managem	ent	127
Using the Display and Control Unit	Creating User F	Profiles	127
Quick Guide: First Weighing	Activating User	rs	130
User Interface (Touch Screen) 32		ofiles	
Using the Touch Screen	Interfaces		133
Activating/Switching Users	USB Port (PC).		134
System Settings (Menu) 3!	PS2 Interfaces	for Bar Code Scanner or Keyboard	135
Leveling the Balance (Q-Level) 30	Interfaces (RS-	232) 25-pin and 9-pin	136
Configuring Calibration/Adjustment	Configuring Se	rial Ports	139
Timer-controlled Actions	Bluetooth® Inte	erface (COM C, Optional)	142
Displaying Device Information 40	Network Interfa	ace (Ethernet)	144
Alibi Memory 4	Data Output		149
Device Parameters44	Data Input		154
Task Management	Updating the Sof	tware	157
Using Applications with the Factory Settings 57	Error and Status	Messages	159
Creating New Tasks (Configuration) 57	GPL License		160
Combining Applications into One Task 60	Care and Mainte	nance	161
Executing Tasks 67	Shipping the Bala	ance	164
Weighing	Disposal		170
Mass Unit Conversion	Specifications		171
Minimum Sample Quantity SQmin	-	ons	
Individual Identifiers 75	Accessories		189
Density Determination	Declaration of Co	onformity	191
Statistics		l Certificate	
Calculation86	Plates and Marki	ngs	197
Averaging92			

Notes on Using this Manual

- ▶ Please read this entire manual carefully and completely before using the device.
- Read the safety precautions carefully.
- ▶ This manual is part of the product. Keep it in a safe and easily accessible location.
- ▶ If the manual should be lost or misplaced, please contact Sartorius for a replacement or download the latest manual from our website: www.sartorius.com

Symbols and Signs

The following symbols are used in this manual:



Warning symbol for various types of dangers.

These symbols are explained in more detail in the section on safety precautions.



This symbol indicates useful information and tips.



This symbol indicates notes on use of the balance in legal metrology within the scope of validity of Council Directive No. 90/384/EEC, replaced by 2009/23/EC (Models MS...-. CE...).



The hand symbol indicates operating steps that involve touch screen controls.



This symbol means that the USER key should be pressed.



This symbol means that the TASK key should be pressed.

Next

When individual buttons are displayed, they should be pressed.

- Indicates a required action
- Describes the result of an action
- 1. If a procedure has multiple steps...
- 2. ... the steps are numbered consecutively.
- Indicates an item in a list



Technical advice / hotline:

Phone: +49 0551.308.4440 Fax: +49 0551.308.4449

Safety Precautions

This device complies with European Council Directives as well as international regulations and standards for electrical equipment, electromagnetic compatibility, and the stipulated safety requirements. Improper use or handling can, however, result in damage and/or injury.

This device should only be operated by trained personnel.

Operators must read these installation instructions, particularly the safety information, and must be familiar with the operation of the equipment. The operator is required to supplement these safety precautions as appropriate. The operating personnel must be provided with the appropriate training.

The requirements pertaining to applicable installation regulations must be followed when using electrical equipment in systems and environmental conditions with increased safety requirements. Relevant laws, standards, regulations, guidelines and environmental protection laws valid in your country must be observed.

Always keep the equipment and balance freely accessible. Any installation work or balance operation that does not conform to the instructions will result in forfeiture of all claims under the manufacturer's warranty.

Danger of Explosion!



Do not use this equipment in hazardous areas in which explosive materials are present.

Danger of Electrical Shock!

If the device is not used in accordance with the instructions, this can affect the product liability for the device.



Make sure that the voltage rating printed on the AC adapter is identical to your local line voltage. The wall outlet must have a protective grounding conductor. The operating instructions included with the AC adapter must be followed.



Do not switch on the equipment if the AC adapter or power cord are damaged.

If the balance itself, the AC adapter or any cables on the balance are damaged, disconnect the equipment from the power supply and secure it so that it cannot be turned on.



Chemicals (e.g. gases or liquids) that can corrode and damage the inside or outside of the balance, AC adapter, power cable or peripherals must be kept away from the equipment.



Do not operate the balance unless the housing and the display are undamaged and the housing is closed so that inadvertent contact with parts inside the equipment is not possible. Make sure that no liquids penetrate the balance and do not allow conductive parts to touch the balance.



Note on Installation:

The operator shall be solely responsible for any modifications to the equipment and for connecting any cables or equipment not supplied by Sartorius.

Information on operational quality is available upon request from Sartorius.

You should only use peripherals and options supplied by Sartorius.

Note on cable quality:

CE conformity is only guaranteed with well shielded cables.

The shielding braid of the cable must be made of metal with at least 80% coverage and also covered with metal foil as much as possible.

The shielding on both sides on the full metal or metalized connector casing, as well as on the casing inlet, must be high-frequency compliant and have a flat design.

You can only turn off power to the device by pulling the plug or disconnecting the power cord, or by unscrewing the four-pin plug at the DC supply cable for the balance.

Do not exert mechanical pressure on the display with sharp objects, as this will damage the display. Liquid penetration can damage the electrical components.

Only use a lightly-moistened cloth for cleaning. See "Care and Maintenance" for cleaning the device. Do not compromise the IP protection of the balance.

Do not open the balance housing. If the seal is broken, this will result in forfeiture of all claims under the manufacturer's warranty.

The device may only be opened by specialized personnel trained by Sartorius.

Disconnect the balance from the wall outlet prior to moving the device.

Avoid condensation damage to the device. When the balance is turned on, condensation build-up is not possible. If the device is moved from a cool environment to a warm environment, you must allow the device to acclimatize for approx. 2 hours before connecting to the power supply.

Hazards at the Place of Installation and During Operation

Avoid generating static electricity and establish equipotential bonding.

A 10 kOhm ground is used when connecting the balance to the power supply. Disconnecting the ground conductor is not permitted.



Exercise caution to avoid glass breakage.

There is a danger of lacerations or cuts caused by sharp edges on glass or metal (contact springs), particularly when removing or replacing the glass draft shield panels; during cleaning; if a glass sample container breaks, or when modifying the balance hardware configuration.



Models with a separate electronics module (e.g. MSx6.6/3.6/2.7):.

Do not replace the electronics module.

Replacing with, e.g. models of the ME/SE series will damage the device.



Make sure all equipment operators have received proper instructions when working with hazardous or toxic materials. Failure to follow correct handling procedures can result in breakage of parts or spillage of liquids or other substances. Make sure the necessary protective clothing or equipment is worn, such as protective gloves, clothing, eyewear. All safety regulations applicable in the laboratory must be observed.



Exercise caution to avoid squeezing or crushing fingers when

- Operating the electronic draft shield
- Removing or installing the panels
- Adjusting the display

Do not expose the device to aggressive chemical vapors or to unnecessarily extreme temperatures, moisture, shocks, or vibration.

The device should be set up in areas away from equipment or devices that have strong magnetic properties. Extreme electromagnetic fields from electrical lines should also be avoided.

Intended Use

Cubis models are high-resolution balances. They were specially developed for exact determination of material mass in liquid, paste, powder or solid form.

Appropriate containers must be used for each type of sample material.

Specific models cover specific weighing ranges: see "Technical Specifications."


Cubis models are designed specially for use in research, education and day-to-day laboratory tasks in science and industry.

They are designed to be used exclusively indoors.

Cubis models can be operated as standalone, connected to a PC, or on a network.

General View of the Equipment and Equipment Supplied

Balances with a readability of ≤ 1 µg

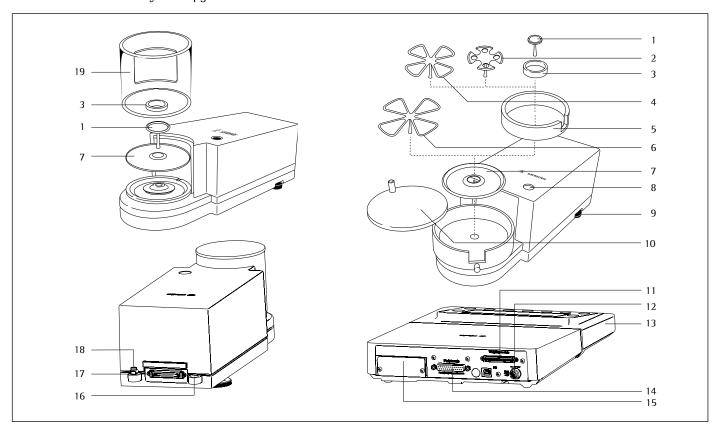


Pos. Description

- 1 Weighing pan
- 2 Filter weighing pan \varnothing 50 mm
- 3 Internal draft shield (for MSx2.7S-F only)
- 4 Optional weighing pan Ø 75 mm
- 5 Shield disk
- 6 Optional weighing pan Ø 90 mm
- 7 Shield plate
- 8 Level indicator
- 9 Leveling foot
- 10 Draft shield cover

Pos. Description

- 11 Female connector for weigh cell
- 12 DC jack
- 13 Display and control unit
- 14 Serial communications port (PERIPHERALS)
- 15 Slot for optional interfaces
 - (e.g., 9-pin data output and PS2 or Bluetooth)
- 16 Lug for attaching antitheft locking device
- 17 Equipotential bonding conductor terminal
- 18 Equipotential bonding conductor terminal
- 19 Draft shield

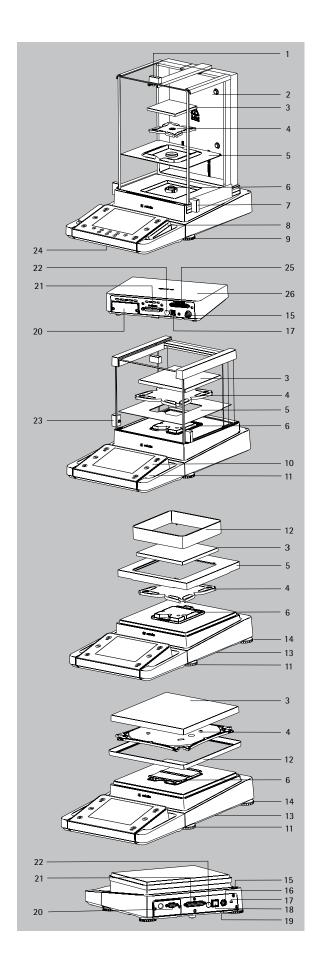

General View of the Equipment and Equipment Supplied

Balances with a readability of ≥0,01 mg

Pos. Description

- 1 Upper sliding draft shield panel/Handle
- 2 Back panel
- 3 Weighing pan
- 4 Pan support (not for MSx225.../MSx125... models)
- 5 Shield plate
- 6 Pan retainer
- 7 Right sliding panel/Handle
- 8 SD card slot (for MSU, MSA models)
- 9 Leveling foot
- 10 Display
- 11 Display and control unit
- Draft shield/Shield disk (only for models with a readability of 1 mg and 10 mg)
- 13 Below-cell weighing port (on the bottom of the balance)
- 14 Leveling foot
- 15 Level indicator
- 16 Power socket
- 17 Slot for attaching an anti-theft device
- 18 USB socket for a PC connection
- 19 Below-cell weighing hook
- 20 Slot for optional interfaces, e.g., 9-pin data output and PS2 (as shown) or Bluetooth
- 21 Communication port (PERIPHERALS) for accessories
- 22 Menu access switch
- 23 Left sliding panel/handle
- 24 Ethernet Interface
 - (on the bottom of the display and control unit)
- 25 Semi-microbalances: Female connector for weigh cell
- 26 Semi-microbalances: Electronics module (for MSx225.../MSx125... models)


Not shown:

- AC adapter
- USB cable
- Operating instructions


Getting Started


Unpacking the Equipment

- Open the box.
- Use both hands to lift the balance, with the packaging, out of the cardboard box.

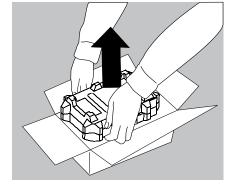

- ▶ Place the packaging with the balance on the floor.
- ► Remove the top part of the packaging.

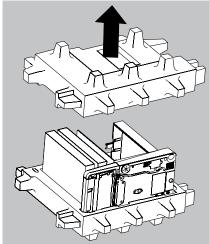

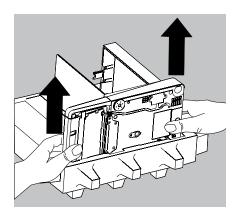



Balance with Draft Shield

- ► Remove the packages (containing draft shield panels, weighing pan, pan support, AC adapter, etc.) from the lower packaging and place them to one side.
- ▶ Use both hands to lift the balance out of the packaging.



Exercise caution to avoid glass breakage. Never lift the equipment by the draft shield panels to remove it from the packaging.

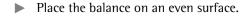


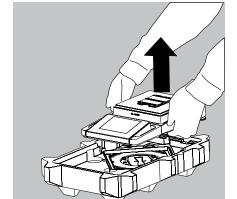

Balance with No Draft Shield

Use both hands to remove the model without a draft shield from the packaging.


▶ Remove the AC adapter and the balance parts from the packaging.

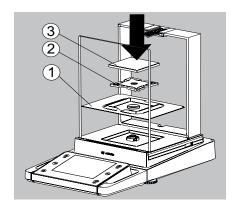


Assembly

Installing the Analytical Draft Shield, Assembling the Balance (Draft shield designation: DA, DI, and DU)


- ▶ Fit parts onto the balance as shown in the picture.
- 1. Shield plate
- 2. Pan support (not for MSx225.../MSx125... models)
- 3. Weighing pan

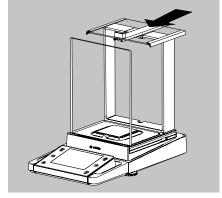

▶ Slide the upper draft shield panel into the guide rails from the rear.

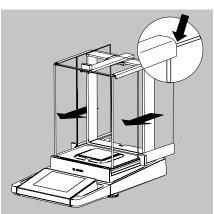


- ► Slide the left and right draft shield panels into the guide rails from the rear. Make sure the panels are within the upper and lower guide rails.
- ▶ Slide the panels in until they engage.
- 1. Right draft shield panel
- 2. Left draft shield panel
- ➤ This completes the balance assembly.

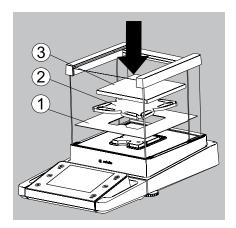


Installing the Draft Shield on Milligram Balances, Balance Assembly (Draft shield designation: DE)

- ▶ Fit parts onto the balance as shown in the picture.
- 1. Shield plate
- 2. Pan support
- 3. Weighing pan

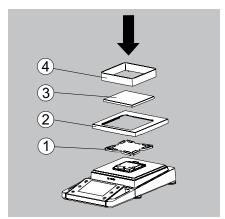

Assembly of Milligram Balances without Glass Draft Shield (DR Option): Balance Assembly


- ► Fit all components listed below onto the balance in the order given:
- 1. Pan support (see also next section)
- 2. Shield plate
- 3. Weighing pan
- **4.** Draft shield frame
- First insert the shield plate.

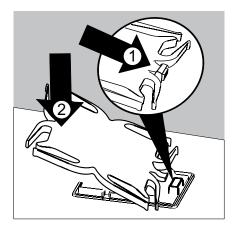


- Make sure the pin on the pan support is facing toward the front, and slide it under the clip on the pan retainer.
- 1. Insert the pan support.
- ▶ After inserting the pan support, press it down against the shield plate so that it lies flat.
- 2. Press down the pan support.




Exercise caution to avoid pinching or crushing fingers.

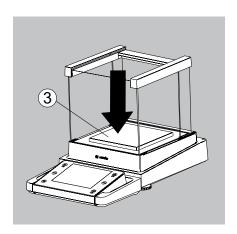


- Place the weighing pan on the pan support.
- **3.** Weighing pan

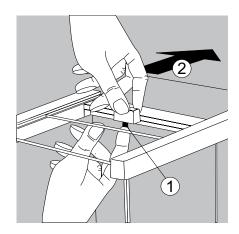

► Slide the upper draft shield panel into the guide rails from the rear while pressing the locking tab.

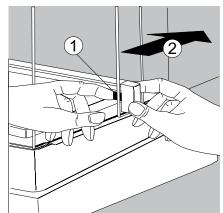


- Slide the side draft shield panel into the guide rails from the rear while pressing the locking tab.
- 1. Right draft shield panel
- **2.** Left draft shield panel

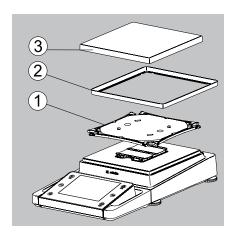

- ➤ This completes the balance assembly.
- ► Remove the panel:
- **1.** Press on the locking tab.
- **2.** Remove the panel
- ▶ Place the panel in the storage slot at the back of the balance.

Assembly for Models without a Draft Shield


- Fit parts onto the balance as shown in the picture.
- 1. Pan support
- 2. Shield plate/Draft shield
- 3. Weighing pan

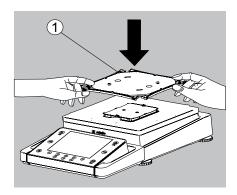

- Place the pan support diagonally and press down lightly.
- 1. Pan support




Carefully turn the pan support clockwise until the two buttons engage. The pan support is now attached.

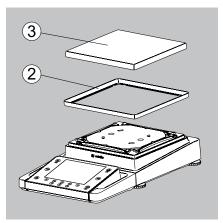

- 2. Insert the shield plate/draft shield.
- ▶ **3.** Place the weighing pan on the pan support.
- ➤ This completes the balance assembly.

MSx225..., MSx125:

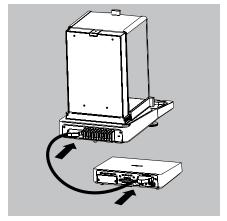


Connecting the Weigh Cell to the Electronics Module/Evaluation Unit

▶ Connect the weigh cell and electronics module using the connection cable.




Please check the plug contacts to ensure a proper connection.

MSx6.6S/3.6P/2.7S:

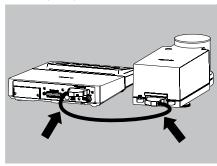

► Connecting the Weigh Cell to the Evaluation Unit

MSx6.6S/3.6P/2.7S:

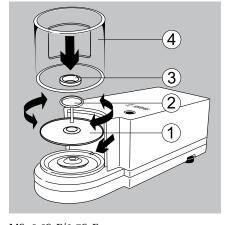

Balances with a readability of 1 μg

Install the following parts in the order listed:


- ▶ 1. Shield plate
- ▶ **2.** Weighing pan

Note:

After inserting the weighing pan, turn it slightly to the left and right, while pressing it down lightly.


- ▶ 3. Internal draft shield (for SE2 only)
- ▶ **4.** Draft shield: Center the hole over the pan (see arrows)


MSx6.6S-F/2.7S-F:

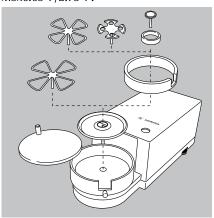

MSx6.6S-F/2.7S-F models

- ▶ Place the components listed below inside the weigh cell in the order given:
- Shield plate
- Internal draft shield
- Filter weighing pan Ø 50 mm or weighing pan
 (optional filter weighing pan Ø 75 mm or Ø 90 mm)
 Note: After inserting the pan, turn it slightly to the left and right, while pressing it down lightly.

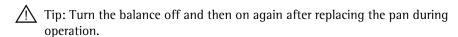


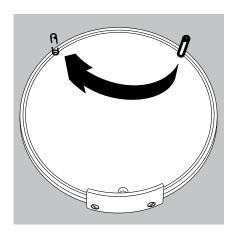

- Internal draft shield (for MSx2.75S-F only)
- Setting up a filter weighing pan for left-handed users:
- Remove the draft shield cover

Draft shield cover


Unscrew the pin and move from the right to the left side

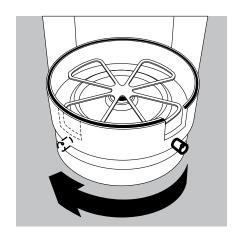

► Turn draft shield parts approx. 90 degrees to the left (loosen knurled screw)


Moving the Balance

Transporting the Device over Short Distances

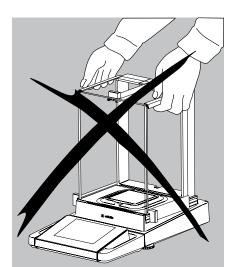


Exercise caution to avoid glass breakage.

Never lift and carry the balance by its draft shield.

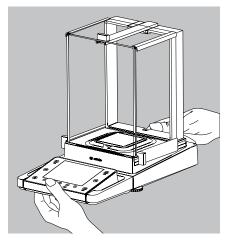

► Carry as shown in the illustration.



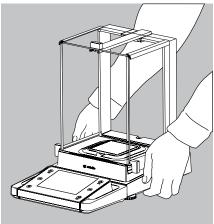

Carry as shown in the illustration.

Storage and Shipping Conditions

- Permitted storage and shipping temperature: -10 to +60 °C
- Unpacked devices can lose their precision if subject to extreme vibrations.
- Excessive vibrations may compromise the safety of the equipment.
 Save the packaging for any future storage or shipment of the balance.
 Only the original packaging provides optimum protection for the equipment.
- Follow all warnings and safety precautions.
- Follow the instructions in the chapter: "Transporting the Balance."

Installation Instructions

- Select the proper setup location:
- Set up the device on a stable, even surface that is not exposed to vibrations (e.g. weighing stone).
- Place the device in a location with enough free space around it so that excessive heat cannot build up.
- Maintain free access to the device at all times.

Choose a location that is not subject to the following negative influences:


- Heat (heater or direct sunlight)
- Drafts from open windows, AC systems and doors
- Extreme vibrations during weighing
- Excessive moisture

Acclimatization

Condensation from humidity can form on the surfaces of a cold device when it is brought into a substantially warmer area.

To avoid the effects of condensation, condition the balance for about two hours at room temperature, leaving it unplugged from the power supply.


Carefully read all warnings and safety precautions in the respective section of this manual.

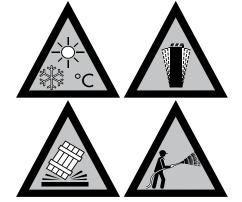

Setting the Angle for the Display and Control Unit


The angle of the display and control unit can be adjusted by the user, in order to ensure optimal readability of the weight values on the display at all times.


▶ On MSA and MSU models, the display and control unit can be tilted as desired.

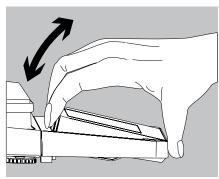

Power Cord Receptacle


- Check the voltage rating and plug design.
- Make sure that the voltage rating printed on the manufacturer's ID label is identical to that of your local line voltage.
- If the stated supply voltage or the plug design of the power cord does not comply with your country's standard, please inform the nearest Sartorius representative or your dealer.
- The power connection must be made in accordance with the regulations applicable in your country.
- In order to connect the equipment to the power supply (protection class 1), use a correctly installed wall outlet with a protective grounding conductor (PE) and a fuse of maximum 16 A.
- If a main feeder from the ceiling is required or if installation of a CEE plug is necessary, these must be carried out by a qualified electrician.
- Only use original Sartorius AC adapters.

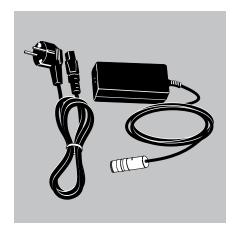

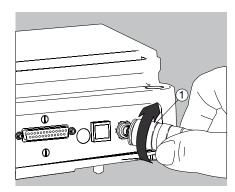


Analytical and precision balances:

1. Plug the DC supply cable of the AC adapter into the power socket of the balance and tighten the threaded fitting.

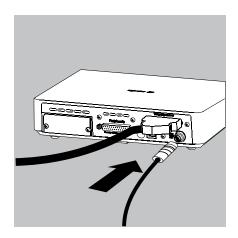

Balances with a readability of ≤ 0.01 mg


1. Plug the DC supply cable of the AC adapter into the power socket of the electronics module and tighten the threaded fitting.

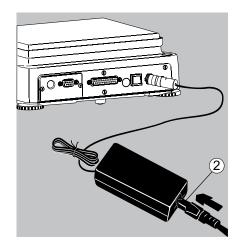

- **2.** Plug the other end of the power cord into the AC adapter.
- **3.** Insert the AC adapter power plug into the wall outlet.
- ➤ The balance is now ready to use.

Safety Precautions

The output line of the adapter has a terminal (GND) connected to the metal housing of the balance. The data port is also galvanically linked to the balance housing (GND).

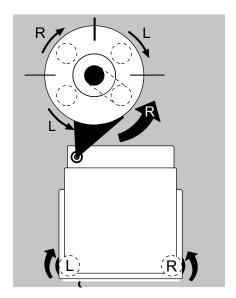

Leveling the Balance, Setting the Level Indicator



Leveling the balance compensates for slant or unevenness at the place of installation. The balance must be perfectly horizontal to ensure consistent, reproducible weighing results. All models are equipped with an electronic tilt angle detection feature. If the balance is not level, all balance models issue an alert on the display.

Push-button automatic leveling is available on models with motorized leveling feet. For models with manual leveling feet, please follow the instructions on the display.

Leveling the Balance Manually

- ► The balance is leveled using both front leveling feet.
- Screw in both back leveling feet (only for models with back leveling feet).
- Turn the two front feet as shown in the illustration until the air bubble is centered within the circle of the level indicator.
- ➤ Normally, several leveling steps are required.
- Screw out both back leveling feet until they touch the setup surface (only for models with back leveling feet).




Warm-up Time

> To deliver precise results, the balance must warm up for at least 30 minutes after initial connection to the power supply. Only after this time will the device have reached the required operating temperature.




When a verified balance used in legal metrology (legal-for-trade applications) is connected to the power, it must warm up for at least two hours before operation.

► Turn on devices via the (♂) key. The following appears in the display:

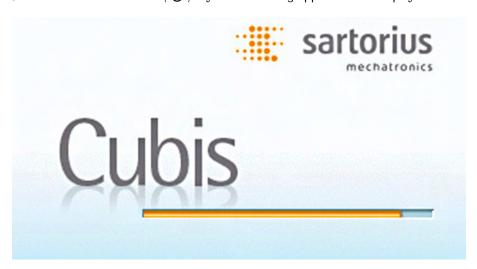

➤ You can now follow the brief instructions to configure balance settings before beginning weighing operations.

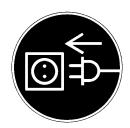


Connecting Electronic Devices (Peripherals)

▶ Make absolutely sure that the device is unplugged from the power supply before connecting/disconnecting any peripheral device (printer, scanner, PC) to or from the data port.




A device connected to the power supply should never be opened.

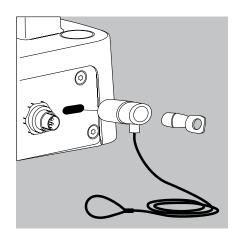

Anti-theft Locking Device (Accessory)


Balances with a readability of ≤0.01 mg

► An anti-theft locking device can be installed to the back of the balance if required.

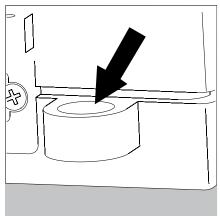

Balances with a readability of $\leq 1 \mu g$


To fasten an antitheft locking device, use the lug located on the back of the weigh cell.

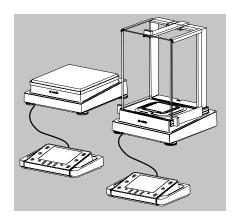

Modifying the Balance


Setting Up the Display and Control Unit at the Place of Use

The display and control unit can be removed for all models to enable the operator to customize the work space.

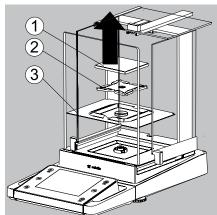

Removing the Retainer with the Display and Control Unit

- Remove all items (such as weights) from the draft shield.
- ► Carefully remove all parts as shown in the illustration.
- 1. Weighing pan
- 2. Pan support (not for MSx225.../MSx125... models)
- **3.** Shield plate/Draft shield
- ► Keep all parts in a safe place.

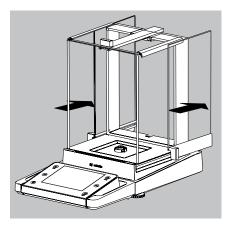


► Remove the panels (right and left).


Models with the analytical draft shield:

- **1.** Press on the locking tab.
- **2.** Remove the panel.


Models with the milligram draft shield:


- **1.** Press on the locking tab.
- 2. Remove the panel.

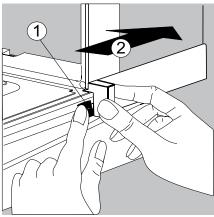

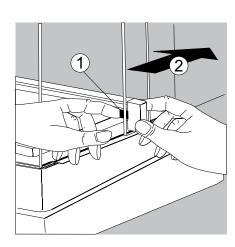

► Keep all parts in a safe place.

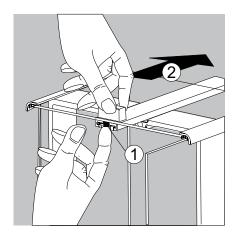

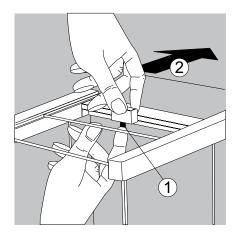


Remove the upper draft shield panel.

Models with the analytical draft shield (left illustration).

- **1.** Press on the locking tab.
- **2.** Remove the panel.


Models with the milligram draft shield (right figure):

- **1.** Press on the locking tab.
- **2.** Remove the panel.
- Keep all parts in a safe place.
- Turn over the balance and place it on a soft surface.

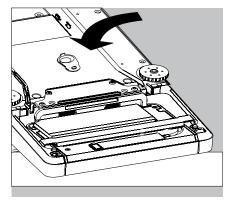





Exercise caution to avoid breaking the glass on models with a draft shield.

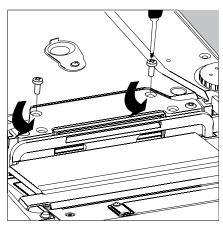

- ▶ Use a 2.5 mm Allen wrench to remove the two screws from the display and control unit retainer bracket.
- ▶ Remove the display and re-insert both screws back into their holes.
- ▶ Lengthen the cable and position the display and control unit as desired.
- ► Turn the balance over and place it on an even surface.

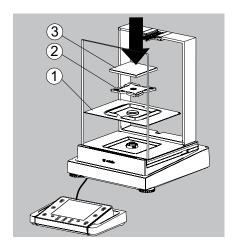

- Carefully place all parts on the balance.
- 1. Shield plate/Draft shield
- 2. Pan support (not for MSx225.../MSx125... models)
- 3. Weighing pan

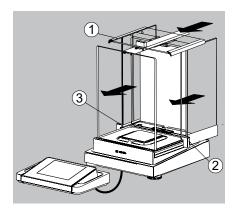


- Replace the upper and side shield panel.
- 1. Upper draft shield panel
- 2. Right draft shield panel
- 3. Left draft shield panel
- ▶ Level the balance.




Leveling the balance compensates for slant or unevenness at the place of installation. The balance must be perfectly horizontal to ensure consistent, reproducible weighing results. All models are equipped with an electronic tilt angle detection feature.

If the balance is not level, the balance issues an alert (MSA and MSU models, see brief instructions). The display shows a graphical representation of a level indicator and text prompts that guide the user through the leveling process.

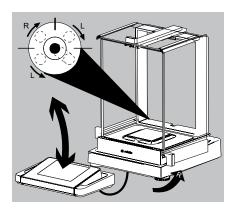


- The balance is leveled using both front leveling feet.
- Screw in both back leveling feet (only for models with back leveling feet).
- Turn the two front feet as shown in the illustration until the air bubble is centered within the circle of the level indicator.
- ➤ Normally, several leveling steps are required.
- Screw out both back leveling feet until they touch the setup surface (only for models with back leveling feet).

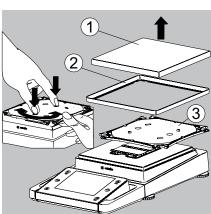


Removing the Display and Control Unit for Models without a Draft Shield


- ► Carefully remove all parts as shown in the illustration.
- 1. Weighing pan
- 2. Shield plate/Draft shield
- 3. Pan support
- ► Keep all parts in a safe place.

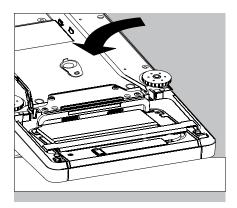


► Turn over the balance and place it on a soft surface.

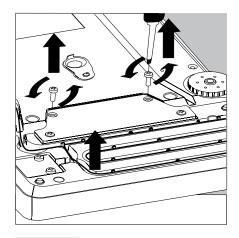


- ► Remove the two retaining screws.
- ▶ Remove the display and re-insert both screws back into their holes.
- ► Carefully pull the cable connected between the display and control unit from the retainer.

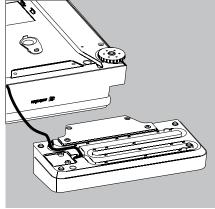

▶ Determine the required cable length.

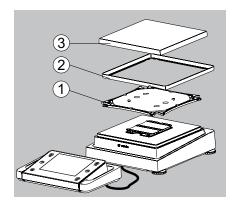

- ▶ Return the balance to an upright position and fit the parts onto the balance.
- 1. Attach the pan support.
- 2. Shield disk (only for models with a readability of 10 mg)
- 3. Weighing pan

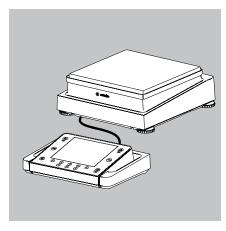

Level the balance.


Semi-microbalances: Attaching the Display and Control Unit to the Electronics Module (MSx225.../MSx125... models)

The display and control unit can also be attached to the electronics module if required for operation.


Turn over the balance and place it on a soft surface.

Remove the connection cable from the cable channel:


▶ Remove two screws from beneath the weigh cell and detach the plate.

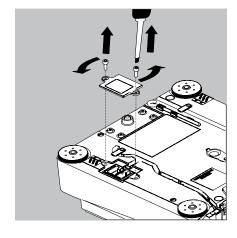


- Remove the connection cable plug.
- ► Then reattach the plate to the slot.

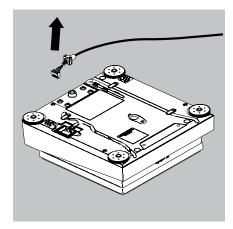


- Remove the display and control unit from the weigh cell: Remove two retaining screws.
- ► Remove the display and control unit.

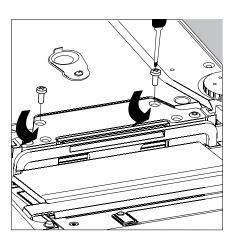

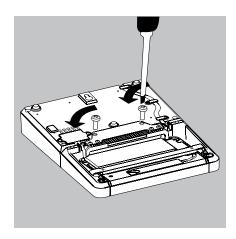


Attach the display and control unit to the electronics module: Reattach the two retaining screws.

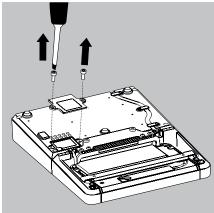


Open the slot for the connection cable on the electronics module:

Remove the screws from beneath the electronics module and detach the plate.

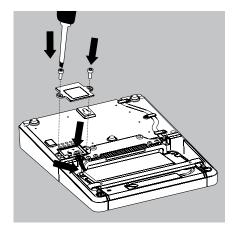

Connect the display and control unit to the electronics module:


- Plug in the connection cable.
- Then reattach the plate to the slot.
- ▶ Slide the cable that sticks out into the cable channel.

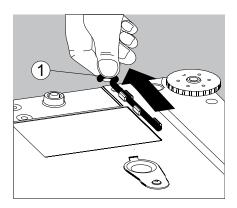


Below-Cell Weighing

A port for a below-cell weighing hook is located on the bottom of the balance.

1. Remove the hook for below-cell weighing from the clip on the bottom of the balance.

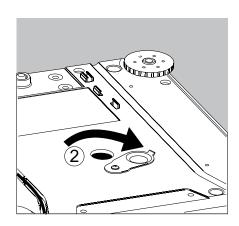

2. Push the cover of the below-cell weighing port to one side.

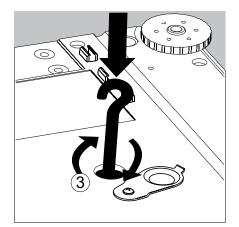

3. Carefully screw on the below-cell weighing hook.




Do not screw it in too tightly, as this could damage the thread or the balance.



Install a draft protection shield.


Attach the sample (e.g., using a suspension wire) to the hook.

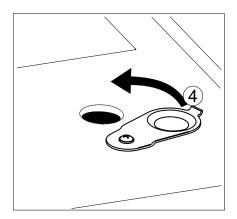

- 4 When weighing has been completed, unscrew the hook and return it to the clip.
- Close the cover of the below-cell weighing port.



The below-cell weighing port may not be opened or used on balances used for legal metrology.

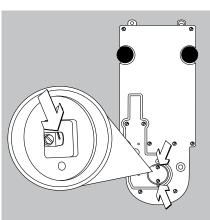


Balances with a readability of ≤ 1 µg

- Remove both screws from beneath the weigh cell and detach the cover plate.
- Attach a wire, for example, to the sample and hang it on the notched hook.




Install a draft protection shield.

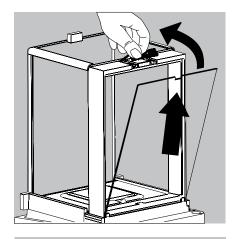


Using the Cable Opening of the Draft Shield

Models with an analytical draft shield have an opening for passing a cable (for example, for a temperature sensor) through to the interior of the weighing chamber.


- 1. Lift the locking tab on the rear panel of the analytical draft shield.
- **2.** Lift the panel out of the draft shield.

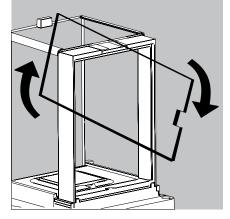

► Turn the panel clockwise 180° so that the opening is at the bottom.

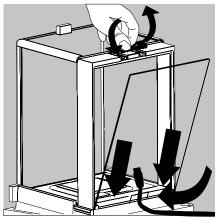

- lnstall the desired sensor.
- Insert the panel in the guide rail.
- ▶ Lift the locking tab and gently press the panel into position.
- ▶ Press the locking tab down and close it.
- ► You can now begin weighing.

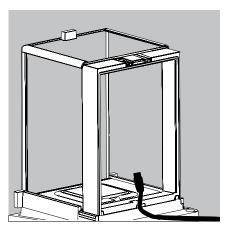

Using the Balance

Turning the Balance On/Off

- ▶ Make sure the balance has been installed and put into operation in accordance with the installation instructions.
- (\bigcirc) Press the on/off key (\bigcirc) on the control unit.
 - ▶ The start screen appears briefly on the display, then the user interface appears.
- On subsequent starts, the most recently active user profile and task are opened (if at least one user profile has been set up).
 - ▶ If prompted, level the balance (see "Leveling the Balance" for details).
- (\circlearrowleft) To put the balance in standby mode or switch it off, press the key (\circlearrowleft).
 - Close the draft shield (if present on your model).

Operating Design: Q-Guide


The Cubis precision and analytical balances are controlled by application software with interactive operator guidance. Once you open a menu or select an application, brief instructions shown on the display will guide you through the menu or application step-by-step. At each step along the way, the display shows only those options that are relevant to your process; this helps prevent unnecessary "detours" and enables you to reach your goal more quickly.

This type of user guidance is mostly intuitive. Although operation of the balance is also mostly intuitive, this manual includes a section with very detailed, step-by-step instructions where you can learn about all of the available options (see "Creating User Profiles" in the chapter "User Management").

Basic Operational Structure

The most basic functions, weighing and taring, can be carried out as soon as the balance is switched on. The application software is divided into three general areas in which individual settings can be configured:

In **Task Management (TASK)**, task profiles with specific applications, weighing and printout settings can be configured. **User Management (USER)** lets you set up user profiles with specific configuration settings and user rights. You can also set up

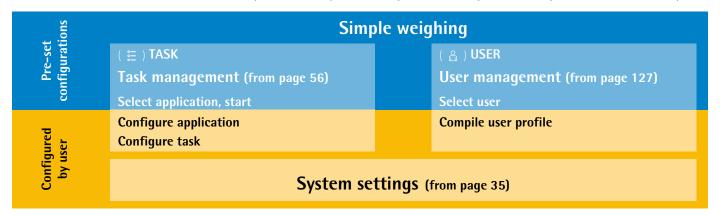

password protection to meet your security requirements. Your balance can also be used for user management.

The **System Settings (Menu)** contains all basic settings and parameters used to operate the balance that affect the entire balance operation. They can also be used for creating new tasks.

The Wizard

You can choose between two display types in several menus:




The overview lists all parameters with their set options. You can select each individual option to change its setting.



When you activate the Wizard in the same menu, the program guides you **step-by-step**: The individual parameters with their options are displayed in succession.

Using the Display and Control Unit

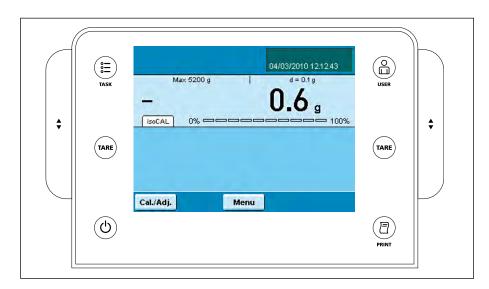

Control keys:




TASK key: Opens Task Management in which applications can be selected and tasks defined. While working in this menu, you can also use the TASK key to cancel at any time and return to task overview.



USER key: Opens User Management in which users can be selected and user settings can be made. While working in this menu, you can also use the USER key to cancel at any time and return to user overview.

TARE) Left and right TARE keys: used to tare the balance.

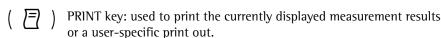

(()) On/Off key: Used to switch on/off and to standby.



Operating elements for electronic draft shield (optional)

Display and Control Unit


The tilt angle of the display and control unit can be set as desired to enable optimal readability for changing working conditions. The color and brightness of the display can also be changed to match the illumination in the room (see "User Management" and "System Settings").

SD Memory Card for Data Exchange


The display and control unit is equipped with a slot for an SD memory card where you can import and export data (task, user profiles). Handling the memory card is explained in chapter **Importing/Exporting Data**.

Setting the Language

➤ The default language for the display is English.

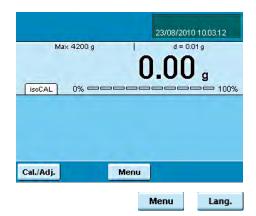

- To change the language, select Menu and then Lang.
- ▶ Use the cursor softkeys to select the desired language.
- > The display changes directly to the selected language.

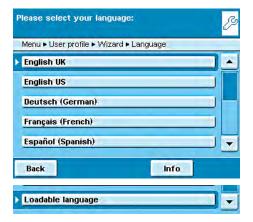

- ▶ If you would like to select a language that was not included with the equipment supplied but was later downloaded from the Cubis website, you must first import the language files (see Section "Importing/Exporting Data").
- To display the downloaded language, select the last option in the language list.


Quick Guide: First Weighing

- (\bigcirc) Press the operating key (\bigcirc) to turn on the balance.
 - Close the draft shield (if available).
- (TARE) Press the tare key (TARE) to tare the balance.
 - ▶ The balance is tared and the displayed value is zero.
 - ▶ If required, open the draft shield and carefully place the sample on to the weighing pan (in a suitable container if required).
 - ► Close the draft shield (if available).
 - ➤ You can read the measured value as soon as the weight value stops changing and the unit is displayed.

User Interface (Touch Screen)

Many operating steps are carried out via the display using the touch screen. The display varies depending on whether the application software is in operating mode or a menu is currently open (setup mode).

Display in Operating Mode

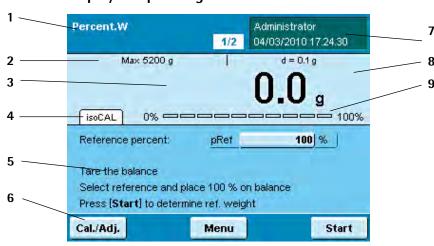



Verified balances feature the following additional displays:

In the line for metrological data:

Min Minimum weight

e Verification scale interval

In the value range:

- o Calculated values.
- pcs Piece count.

- Function display: Current task (here: Step 1 of 2) Touching this area displays the description of the task.
- **Line for Metrological Data:** left: Weighing Capacity (max); right: **d** = Readability (resolution) of the balance.
- Value range with current measurement value (the numeric value is displayed in black numbers once the weighing system reaches stability; beforehand it is displayed in gray numerals)
- Field for status and warning messages (see below)
- 5 Task area with instructions on how to proceed
- 6 Toolbar with currently available buttons
- User field: Displays the current user, date and time Touching this area displays the description of the user.
- For _.CE model balances verified for use as legal measuring instruments: Displays the weights that are not weight values, e. g. calculated values
- Scaled measurement value display (percentage of used weighing capacity)

Status and Warning Messages



tabs with the following messages may be shown here:

isoCAL is displayed: isoCAL is active (extended temperature range)

isoCAL flashes: Calibration is required Touch the tab to activate the function.



Level me flashes: Balance is not level, leveling is required

Touch the tab to activate the function.

GLP: GLP print is active, the header will be printed. Touch the tab to print the footer.



SQmin: The SQmin value entered is permanently displayed.

SQmin flashes: the weight value is less than the entered SQmin value.

U 12.00 g


U or **U*** or **PA**: The selected DKD value (measurement uncertainty) is currently displayed.



Ion: lonizer in progress (only when ionizer available)

Touch the tab to deactivate the function.

Display in Setup Mode

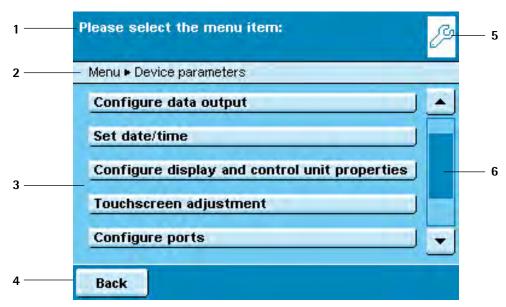

- 1 Interaction area with instructions on how to proceed
- **2** The orientation line shows the current path
- **3** Working environment with selectable options
- 4 Toolbar with the available buttons
- 5 Symbol indicating which menu (TASK, USER or MENU) is currently open
- **6** Scroll bar to page through the selection area

Using the Touch Screen




Equipment may be damaged by objects with points or sharp edges (e. g. pens).

The touch screen should only be operated using the finger tips.




All **operating elements** on the touch screen are depicted in a light color with a shadowed 3D effect. They are activated with a light touch.

Alternative selection: When only one of several options is available, always touch the desired option and the program will automatically switch to the next view.

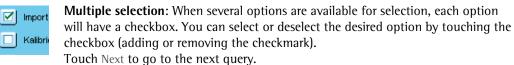



For some displays, you will need to use the **scroll bar** to view all options. You can either slide the bar (touch it and pull in the desired direction) or touch on the up or down arrow keys.

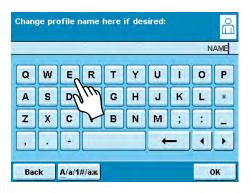

Entering Text and Numbers

A keyboard will appear whenever you have to enter text and numbers. A cursor appears in the entry line above the keyboard.

- Select the desired character with a light touch.
- ➤ The entered text will appear in the entry line.



Press the **shift button** to change the input mode between uppercase letters, lowercase letters, numbers and the complete character set with special characters.



The **backspace key** is used to delete the character to the left of the cursor.



Toggle input mode

Touching the entry line toggles the input mode.

Overwrite mode: The cursor marks an individual character or the entire entry. The marked text will be overwritten by the new entry.

Input mode: The cursor is positioned in the blank space between characters so that you can enter more characters.



Use either of the **Arrow keys** to move the cursor one space to the left or to the right.

Back

The **Back** key cancels the entry process and takes you back one step to the last view

ок

The **OK** key ends the process and saves character input.

Activating/Switching Users

When the balance is in the operating mode, the active user is shown at the top right of the display in the user field. To activate another user, open the User Management menu.



- ► Press the USER key (🖺).
- > The user selection is displayed. The current, active user can be recognized by the small arrow and the dark background.
- Touch the user you want to activate.
- ▶ If password protection activated, enter your password.
- ➤ The program switches back to operating mode, with the selected user profile activated.
- ▶ When you touch Start, the user who is currently selected will be activated.



System Settings (Menu)



This menu summarizes all the balance settings relevant to the device. Basic settings can be made here that immediately affect the device. Changed settings do not affect previously defined tasks and user profiles.




All system settings are user-independent, i. e. they are valid for all users.




As long as a task is active and/or a user has been selected, several basic settings cannot be changed. In order to change all settings, "Weighing" has to be selected in the operating mode and no user may be selected.



Not all functions/settings can be selected on balances used as legal measuring instruments!

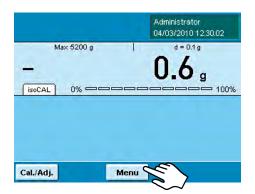

To open System settings, touch **Menu** when the balance is in the operating mode.

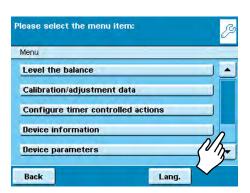

- > The list of available system settings is displayed.
- To view all menu items, use the right scroll bar.
- Touch the menu item whose settings you want to change.

System settings are divided into the following menus:

- Leveling the balance
- Configuring calibration/adjustment
- Timer-controlled actions
- Display device information
- Configure device
- Import/export data
- Activate service mode

You have already learned how to use the menus (see section "Using the Touch Screen"). Only the configuration options and their meanings are described below.

Leveling the Balance (Q-Level)

It is essential for exact weighing results that the balance is absolutely level. The front leveling feet can be used to level out small tilts of the floor. Depending on the model, the balance is equipped either with leveling feet that can be operated manually or motorized leveling feet.

All Cubis balances are equipped with the automatic **Q-Level** feature. An integrated sensor detects the alignment of the balance and triggers a warning message when leveling is required. You can configure the leveling notice and the automatic leveling function (see System Settings/Device Parameters).

When the isoCAL function is activated, the isoCal trigger is set after manual leveling. Balances with motorized leveling feet are leveled automatically prior to each new calibration/adjustment.

Level me

When "Level me" appears on the display, you must level the balance. The process is quick and easy.

Level me


Touch the tab Levelme

٥r

Menu
Level the balance


- Select the option Leveling the balance from the menu.
- Make sure that the weighing pan is empty.

Balance with motorized leveling feet To start the leveling function, touch **Start**. Follow the instructions on the display.

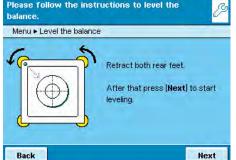



Follow the instructions on the display.

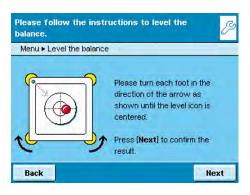



After each leveling routine, the balance must be **calibrated/adjusted** (see chapter **Calibration and Adjustment**).

Configuring Calibration/Adjustment



Menu

Setting options are divided into the following sections:

- Defining calibration/adjustment functions
- Defining external check weights
- Setting isoCAL fully automatic adjustment function
- Enter the desired settings in all submenus.
- To save settings, touch Save.

Factory settings are marked with an *.

Defining Calibration/Adjustment Functions



cal./adj.key: Function of the key: Block key/*Configure key selection menu/ Select fixed 1-key function.

cal./adj. selection menu: Which calibration/adjustment functions should be offered when touching the CAL key?

Cal./adj.sequence: Options: *Calibration, then auto.perform adjustment/calibration and then manual adjustment.

Cal./adj. unit: Options: *gram/kilogram/defined by data record.

Calibration/adjustment data Define external calibration weights

Defining External Check Weights

Number of ext.check weights: Enter the number.

External check weight n: Enter the data for each check weight: Weight value

(in grams)/ID/Certificate/Validation

Calibration/adjustment data Set fully autom. adjustment function isoCAL

Configuring the Fully Automatic Adjustment Function isoCAL

isoCAL: Defining the function: Off/Only warning level, manual trigger/*On, automatic implementation

isoCAL warning level: *Only isoCAL status field/Warning message, repeat/ Alarm message, adjustment mandatory.

isoCAL temperature: Enter temperature differential (in Kelvin).

isoCAL interval: Enter interval in hours.

Timer-controlled Actions

You can program the following actions so that they are run automatically at prespecified times:


- Display a message and execute a command (see example below)
- Execute a command
- Show a message (e. g. a prompt as a reminder)
- The available actions are displayed. When you open this menu for the first time, the list will be empty.
- ► Touch **Edit**.

Menu

Edit


To define a new timer-controlled action, touch **Create**.

Here you can later also modify, copy or delete each action.

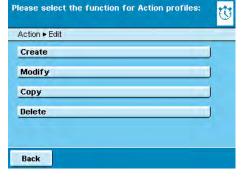


Timer controlled action


Example: You want to have the "Internal cal./adjustment" function performed each morning at 7:00 a.m. In addition, you want have a message displayed.

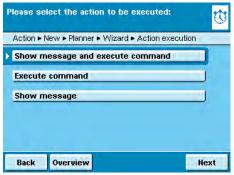


 Select the action you want to install (e. g. Display message and execute command).

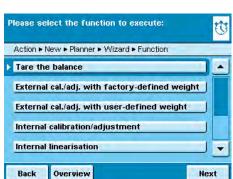

Select the command you want to have executed (e. g. Internal cal./adjustment).

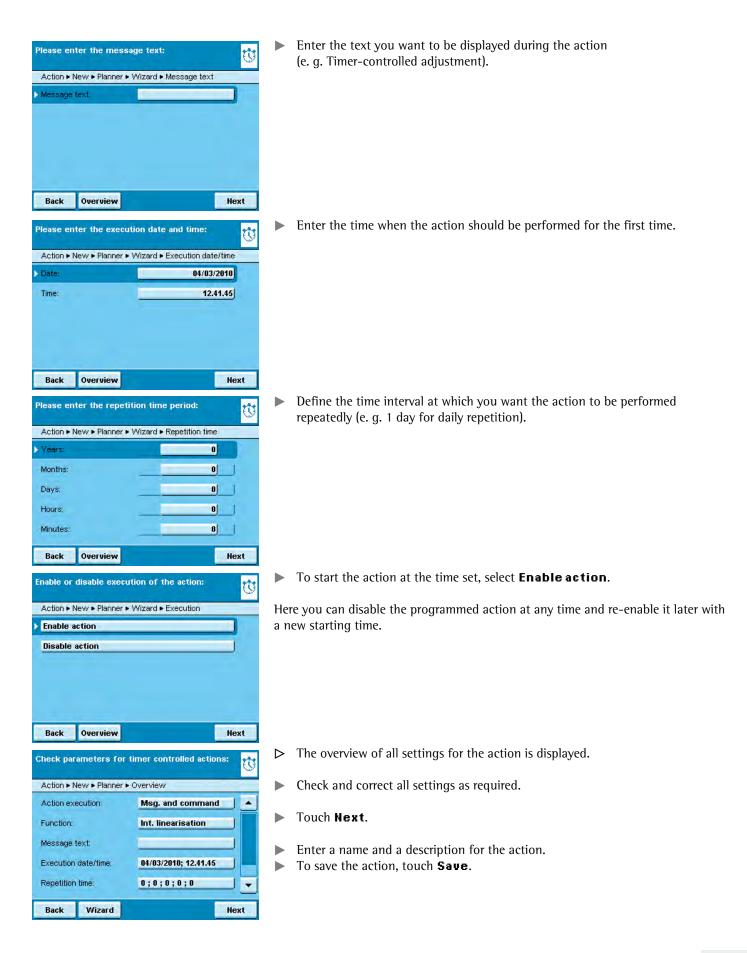

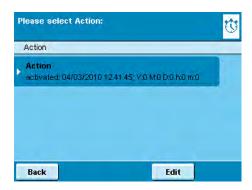

▶ The programmed action is displayed in the list of actions. Here you can edit them again at any time.

Displaying Device Information



You can view various information here.

Select the information that you want to view.

Viewing Basic Information

Here you can find all information about manufacturer, model, serial number, host name and IP address used.


SQminDKD ► To display SQmin and DKD parameters, touch **SQminDKD**.

► To view the version of the balance, display and control unit and application software, press **Version**.

To view the weighing ranges and digits (resolution) touch **Ranges**.

Information about Service and the Hotline

Service contact information is displayed here along with the hotline number as well as maintenance contract and maintenance interval information.

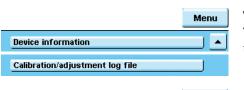

Device information

Service information/hotline

Geräteinformation anzeigen

Audit Trail

Version

Ranges

Menu

Menü

•

Viewing Calibration and Adjustment Data

The log file is displayed here with the results of all calibration and adjustment routines.

Viewing Recordings in Audit Trail In Audit Trail, all changes are logged that have been performed in the Menu, in the user profiles and in the task profiles. The following will be recorded: User, profile, parameter ID, parameter, old and new setting.

Alibi Memory

The alibi memory contains (stores) weighing results with date and time and process number. This takes place every time a print event is triggered (e. g. by pressing $\mathbf{Print}(\ \ \ \ \)$). Tare and input values are also saved. The alibi memory can therefore be used as a $\mathbf{replacement}$ for a \mathbf{log} printer.

The data memory has a ring buffer. The oldest data records can thus be overwritten by new ones. The user can define a memory-time dependent overwrite protection as well as a storage period if desired (default: 90 days, see Section "Configure Device"). The alibi memory can contain approximately 450,000 data records, the storage capacity depends on the volume of the individual data records.



The alibi memory is verifiable. When the balance is verified, you must check the correct function of the alibi memory by saving some values.



If the alibi memory is used in legal metrology (according to Council Directive 90/384/ EEC, replaced by 2009/23/EC, you should observe the following instructions:

- Make sure that there is sufficient storage capacity. Responsibility for this lies with the operator.
- The weighing results should be stored together with a unique ID code (e. g. number or date and time). This makes it possible for every weighing operation or weighing result to be assigned and verified if necessary without any problem. These ID codes must be indicated on any receipts generated on peripheral devices.
- When data are output, all the required tare and net values will be saved.
- It must be possible for the contractual partners involved to be able to check and verify the stored weighing results.
- Any continuously indicated weighing results that are not previously stored and have no identification codes may only be used in applications that are not for use in legal metrology (e.g. batching, controlling or the like). It is not permissible to circulate any results lacking identification codes for applications used in legal metrology!
- Receipts can be prepared by a peripheral device not subject to legal verification according to the preamble to Annex I of Directive 90/384/EEC (2009/23/EC) and Section 7b (3) 2 of the German Verification Act.

These receipts must contain the following information:


- An ID for every weighing result;
- when used in Germany, an additional indication that the weighing results may be compared with the weighing results legally stored on basis of the identification.

You must first install and activate the alibi memory in order to use it (see Device Parameters/Configuring the Alibi Memory).




Searching and Displaying Data in the Alibi Memory

You can search for data using various search criteria and view the storage capacity of the alibi memory.

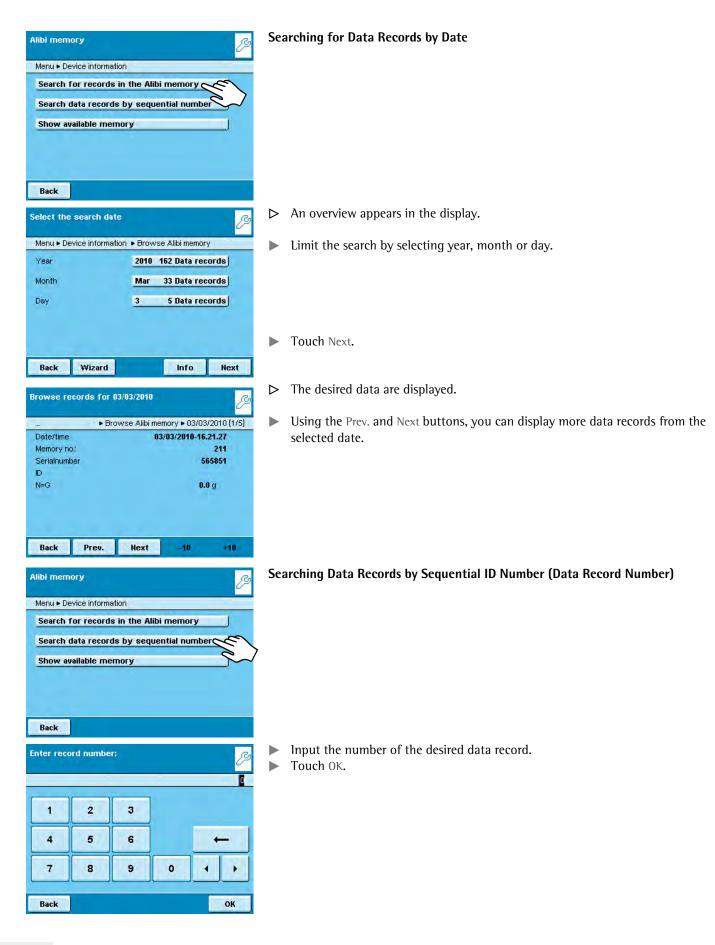




➤ The desired data record appears in the display.

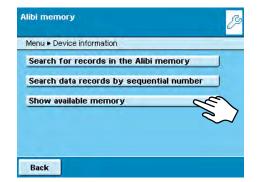

Viewing Information on the Alibi Memory and Available Memory Space




All important status information on the alibi memory are shown here. You can view the percentage of storage capacity utilized and an estimation of the number of data records that can still be stored. The exact number depends on the volume of the individual data records.

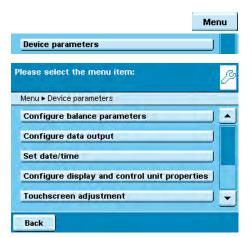

Device Parameters

Here you can change the following equipment settings and update the software. All settings are user-independent.

- Balance parameters
- Data output
- Date and time
- Display and control unit properties
- Touch screen adjustment
- Interfaces
- Alibi memory
- Loading a software update
- Motorized draft shield (optional)
- lonizer (optional)

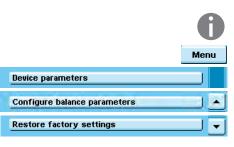

Configuring Balance Parameters

Factory settings are marked with an *.

General Weighing Settings


Here you can configure the standard settings for the general weighing functions. These settings can be changed for a specific task whenever a new task is created.



These settings can only be changed when the balance is set to "Weighing" in the operating mode. If **General weighing settings** cannot be selected, you can touch **TASK** (;) to toggle to Task Management and select **Weighing**.



In order to receive detailed information on the individual setting options, touch **Info**.

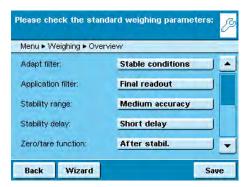

Adapt filter: This setting defines the measuring time of the balance to filter out the effects of unfavorable weighing conditions such as drafts or vibration. Options: Very stable/*stable/unstable/very unstable conditions

Application filter: Equalizing the load deviations in the display. Options: *Final readout/Filling mode/Low filtering/W/o filtering

Stability: The stability indicator lights up when the weight results are constant within a defined range. Options: Maximum precision/very precise/*precise/fast/very fast/maximum speed.

Stability delay: This setting lets you compensate for slowly abating disturbances, such as turbulence in the weighing chamber of an analytical balance). Options: *Very short/short/medium/long delay.

Zeroing/tare function: Conditions for carrying out the tare function. Options: Without/*after/at stability.

Auto zeroing: When this option is turned on, changes of a set fraction of scale intervals per second from the display zero point are tared automatically. Options: On/*Off

Base unit: Weight unit to be used for this task. Options: All available weight units. *gram.

Display accuracy: A lower display accuracy results in a faster display.

Options: *All places on/Last place off

Tare/zero at Power: Automatic taring when the balance is switched on.

Options: *On/Off



Setting up Q-Level Leveling Function

You can configure the leveling function here.

Leveling notice

Save

Off, no notice: The prompt to perform leveling is omitted.

Status display: As soon as leveling is required, the tag **Level me** appears and flashes.

Alarm message, leveling mandatory: As soon as leveling is required, the tag **Level me** appears and flashes. If this notice is ignored, an alarm message appears after 5 minutes. In that case, you can only continue working after you have performed leveling.

Automatic leveling (only on balances with motorized leveling feet.)

Off, start by touching: To activate automatic leveling, touch Start.

Activate the fully automatic function: Before every calibration/adjustment routing, the balance will be leveled automatically.

► To save settings, touch Save.



Back

Defining the Startup Mode of the Balance

You can select the mode that is used when the balance is turned on and when you touch the (🖒) operating key repeatedly.

Options: On/standby or Always on

To save settings, touch **Save**.




Defining the Power saving Mode of the Balance

Switch the power saving mode of the balance on or off. **Off, no function**: The power saving mode is not in use.

Auto shutoff after 2 minutes: The backlighting of the display and control unit switches off if the device is not operated for 2 minutes.

► To save settings, touch **Save**.

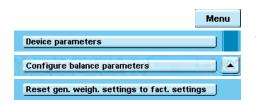

Resetting General Weighing Settings to Default

You can reset all balance settings back to the default delivery status (after a security prompt).

Configuring Data Output (Print Parameters)



Here you can configure all settings for printing and data output. Several of these settings depend on the interface configuration (see chapter Interfaces).

Factory settings are marked with an *.

In order to receive detailed information on the individual setting options, touch **Info.**

Interface to the print output: Options: COM port A, B, C (file/SD memory card), D protocol: Options: Print/SBI/XBPI/Web service/SICS/Remote display. This setting cannot be changed here, but is defined when configuring the interfaces

0

Which parameters and settings are displayed here depends on the configured protocol.

GLP standard print: When should the ISO/GLP output be possible? Options: *Off/enabled for cal.-adjustment/always enabled.

Taring after printout: Automatic tare after printing? Options: *Off/on Print event: What event should trigger printing?

Options: *PRINT key/Start of task/Initialization of application/Result of application/Printing component/Evaluation of application/Exit task/Calibration - adjustment Print key elements: Which elements should be printed?



For further information, please refer to the chapter **Interfaces**, in the section **Configuring a Printout**.


Setting the Date and Time

Device parameters

Set date/time

► Enter the current date and time in the prescribed format.


Configuring Display and Control Unit Properties

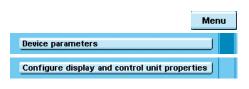

The following properties of the display and control unit can be changed, as necessary: Language, background colors, brightness, acoustic signal.

These general user-independent properties are used when no user is logged in and serve as a template when a new user is created.


Adjusting the TouchScreen

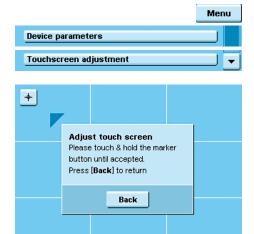

Here you can customize the marker button on the touch screen (finger pressure, angle of touch, etc.).

► Follow the instructions on the display.

Configuring Interfaces and Network Settings

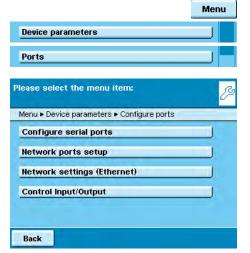


Here you can configure interfaces (serial ports and network) and make settings for the network.

A detailed description of the connection options and data interface ports can be found in the chapter **Interfaces**.

Configuring Serial Ports


Menu

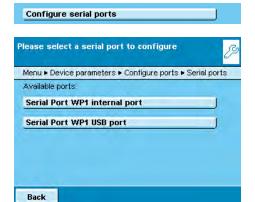

Device parameters

Configure ports

- > The available interfaces are displayed.
- Select the interface you want to configure.
- ▶ Determine all settings for this interface.
- To save settings, touch **Save**.

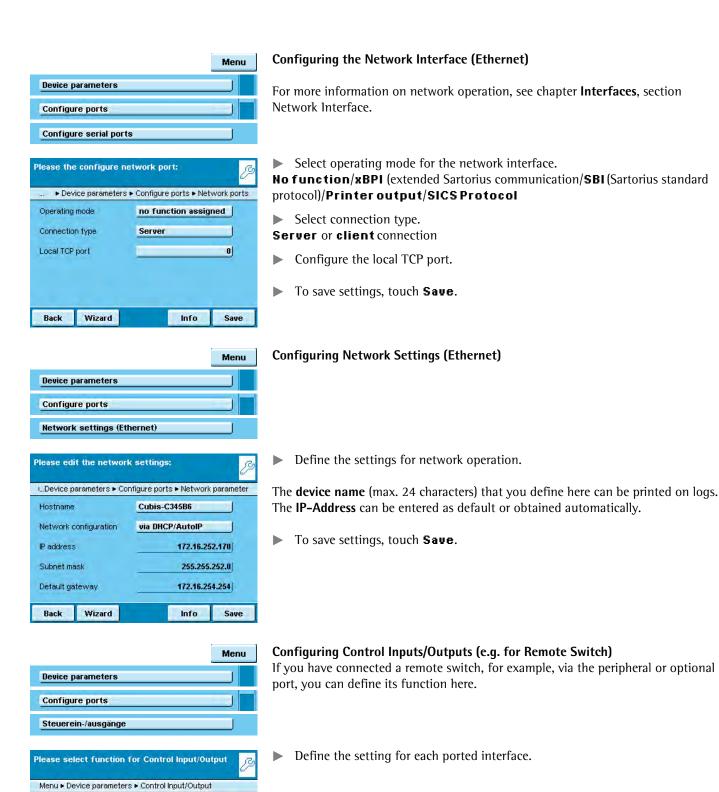


Back


Control inputs/outputs:

Foot switch YPE01RC


Save

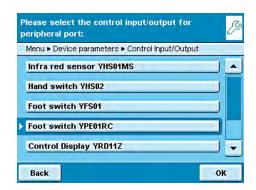

- ▶ Select the ported hardware device or control function.
- ① During checkweighing, the control outputs have no function when set to:
 - Foot switch YPE01RC
 - Control inputs

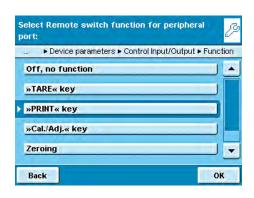


- Define the settings for the control inputs/outputs. Off, no function/TARE key/PRINT key/Cal./adj. key/Zeroing/taring
- ➤ To save the settings, touch Save.

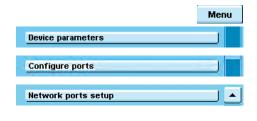


Configuring "Print to file"

Purpose:

- Transferring a print file in TXT or CSV format to a SD card or FTP server.
- Viewing a print file directly in an Internet Browser.

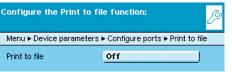

Select Off, no function.

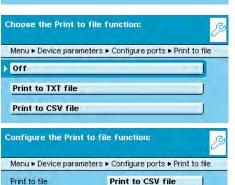

Off

Manual data transfer

Automatic data transfer

Specify the file format.

Print to CSV file:

In addition to the SBI output (like in the TXT file), three additional columns are also written to the file: ID, Value and Unit. The 4 columns are separated by the ";" character. The "+" sign has also been removed in the third column and the decimal point has been replaced by a decimal comma (because of Excel).

- ► Touch **Off**, **no function** to specify the settings for automatic or manual data transfer.
- ➤ The data are deleted from the balance after the automatic or manual transfer so that only the new data are transmitted the next time there is an automatic transfer.
- ▶ Manual transfer:

Touch **Trans.SD** or **Trans.FTP** during operation to transfer data to the SD card in the balance or to the FTP server.


- ➤ The transfer takes place in the data/print directory.
- Select the output medium for the file.
- ➤ This file can be transferred manually or automatically to the SD card of the FTP server. Combinations are also possible, e.g. manual transfer to the SD card and automatic transfer to the FTP server.


"To SD card:"


"To FTP server:"

▶ When exporting, print files can also be exported to the SD card independent of the transfer. This will create a copy (backup) of the file to the SD card. This does not delete the file from the balance. After the export, the files are located in the directory data/all/EXPORTxxxx/print.


- ▶ If you use the FTP server option for the transfer, the parameters for the FTP server must be entered in the menu (see next section). The user must have the rights to create subdirectories on the FTP server.
- ► Select the desired interval, the date and the time of the first transfer.
- To save settings, touch Save.

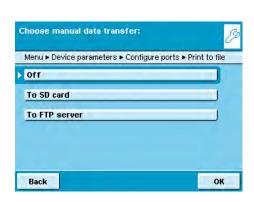

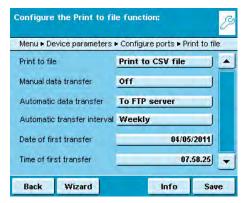

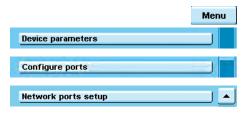

Configuring data transfer settings for a FTP server

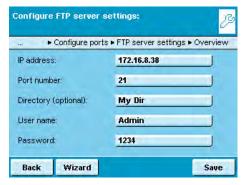

- Specify the data transfer settings to a FTP server.
- Enter the following parameters for the FTP server: IP address, port, FTP user name, FTP password, subdirectory also optional. The user must have the rights to create directories on the FTP server
- For more information about displaying a print record via an Internet browser, go to Interfaces, "Remote Control of the Balance from a PC via Ethernet" section.

Menu Device parameters Alibi memory configuration

Configuring the Alibi Memory

Here you can switch on the alibi memory, delete the recorded data and set the interval for automatic deleting.

Slide the right scroll bar in the submenu **Device parameters** downward to the menu item Alibi config.

Turning the Alibi Memory On/Off Please select configuration Menu ➤ Device parameters ➤ Alibi config.

Alibi memory on/off

Back

Automatic deletion of older data

Reset and initialise Alibi memory

Select whether you want to have the alibi memory switched on or off.



- Define whether the data stored in the alibi memory should be deleted automatically at regular intervals after a pre-specified time.
- As appropriate, enter the desired interval (in days).

Initializing and Switching on the Alibi Memory


You can delete the Alibi memory immediately. The settings for the alibi memory (On/Off, automatic deletion) will remain stored.

To delete data in the alibi memory, confirm the security prompt.

Loading a Software Update

Cubis software is updated at regular intervals. A detailed description of how to perform a software update on your balance can be found in chapter Updating the Software.

Cubis MSA User Manual

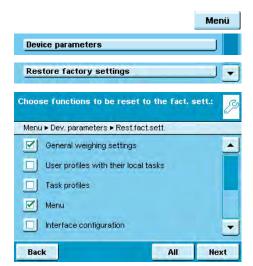

Restoring Factory Settings

You can reset all settings individually to the factory settings at any time.


- Use the cursor softkeys to navigate through the Device parameters submenu to Restore factory settings.
- ▶ Mark any settings you want to reset and select **Next**.
- ► Confirm the reset when prompted.

Installing the Motorized Draft Shield (Optional)

On balances equipped with a motorized draft shield, the control unit has buttons on the right and left sides. These buttons can be used for motor-driven opening and closing of the draft shield doors. You can configure the functionality of these keys in the menu.

Models with the analytical draft shield:

For example, you can configure the left key to open and close the right and upper door (for right handers) and the right key to open and close the left door (for left handers). The opening width (entirely or partially) of the sliding panels can also be defined. You can configure these settings manually on the draft shield itself (see "Learning Function" below).

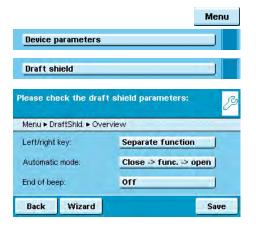


- Check and correct all settings as required.
- To save settings, touch Save.

▶ Define the functions for the left and right keys: Off: The keys have no function.


Same function: Both keys have identical functions.


*Separate functions: The left and right keys are configure with different functions.

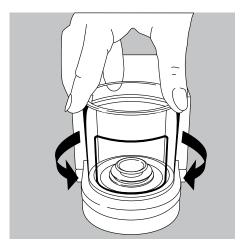

Models with a rotation draft shield (MSx6.6S/3.6P/2.7S):

	Same function (one opening angle)	Separate function (2 opening angles)
Defining the opening width	Manually select the desired opening angle	-
Touch the \$ key	 Open to the previously selected opening angle Close 	1) Touch the left \$\display\$ key: Opens 260° clockwise Right \$\display\$ key: Opens 100° counter clockwise 2) Close

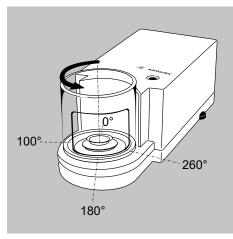

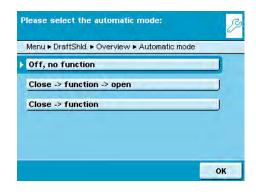

Configure the automatic mode for the draft shield.

*Off, no function: Automatic mode is off.

SClose > **function** > **open**: The draft shield doors are closed in automatic mode while the function is executed, and then the doors re-open.

Close > function: The draft shield doors are closed in automatic mode while the function is executed.

Automatic mode can be executed using the following keys:

- TARE key
- PRINT key
- Cal./Adj. key
- or can be executed when a component is saved for logging applications (statistics, formulation, totalization)



If automatic mode is on, the draft shield doors automatically close when the balance is switched to standby.

Assigning Opening Functions to Keys (Learning Function)

Example: The keys should be cross-configured (left key opens/closes right door and vice versa).

The balance must be turned on.

- Close all draft shield doors.
- Manually operate the door you want to assign to the left key: Push the right door open to the extent you want it to be opened (entirely or partially).
- \triangleright Press the left key \diamondsuit .
- ➤ The door will close again. Now you have configured the left key so that, when pressed, it opens or closes the right door.
- ▶ Proceed accordingly to configure the left door/right key.



According to the same pattern, you can assign operation of the upper door to one of the keys. You can also configure one key to move both doors by manually opening both doors during set up before you press the desired key.

Installing the Ionizer (Optional)



If your balance is equipped with an ionizer module, you can configure the basic settings for the ionizer module here. During any **Electrostatic Charge Eliminator** task, you can change these settings as needed.



Define ionizer function.

Switch off ionizer: The ionizer is switched off.


Manual activation: lonization can be turned on at any time by pressing the **Ion** key.

*Automatic execution after closing draft shield: lonization is executed automatically, as soon as the draft shield is closed.

Installing the Ionizer (Optional)



If your balance is equipped with an ionizer module, you can configure the basic settings for the ionizer module here. During any **Electrostatic Charge Eliminator** task, you can change these settings as needed.

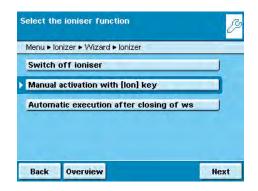


Define ionizer function.


Switch of figures: The jonizer is switched off.

Manual activation: lonization can be turned on at any time by pressing the lon key


*Automatic execution after closing draft shield: lonization is executed automatically, as soon as the draft shield is closed.

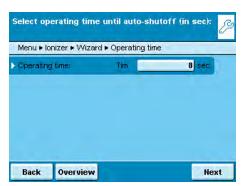

▶ Define how long an ionization routine should take.
Operating time in seconds: 3 to 15 sec, *8 sec.

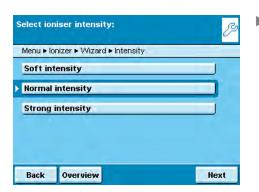

Define the intensity level at which the ionizer operates.
 Low, *Medium, High

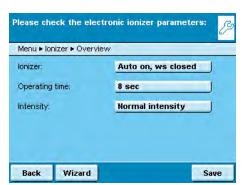

- ► Check and correct all settings as required.
- ► To save settings, touch **Save**.

Task Management



In the Task management menu (TASK), you can:

- Start a task
- Configure a new task
- Edit an existing task (modify, copy, delete)

For **task** configuration, there are numerous **applications** available, e. g. weighing, parts counting, calculation. Each task contains at least one of these applications. However, it can also have a combination of several applications. Each application can be used either with their factory settings or you can also modify them as desired. The **Weighing** application is available as a basic function for every task. Each task also contains the individual balance settings and the individually configured printout.

The options available to you in Task Management depend upon your user access rights: An **Administrator** can use all options while a **User** can only use limited options (see "User Management" in the User Management chapter). If no users have been defined, everything is accessible, there are no rights.

The administrator can **create tasks with restricted use**, i.e. tasks that can only be used by certain users or only by the administrator himself.

Global tasks can be used by all users.

Separate tasks are only accessible to the users who created them.

Exception: A user can copy global tasks to use them afterwards as a local task.

Option	Administrator	User
Configure, modify, delete separate tasks		\square
Configure, modify, delete global tasks	\square	
Execute or copy separate tasks		\square
Execute or copy global tasks	lacksquare	\square







Initially, this overview is empty. Once tasks have been configured, they will be displayed here.

When you touch **Start**, the task currently selected will be started (indentifiable by the dark background).



Importing tasks:

You can also load preconfigured tasks to the balance:


- Unpack the desired ZIP archive file to the root directory (Root Dir) of a SD card after download from the Sartorius homepage or from the included CD.
- 2) In the Balance menu, submenu "Import/Export data": Execute "Import files from the SD card."
- 3) Insert the SD card into the balance.
- 4) Select the corresponding import file and active the task download via **Next**. Then exit the menu.

Using Applications with the Factory Settings

Each application comes with specific factory settings.

If you want to apply these settings unchanged, you can do this quickly during configuration:


Open Task Management and touch Edit.

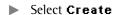





Please select the function for Task profiles:

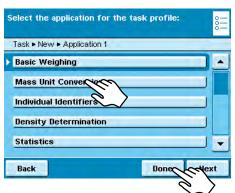

- An application list is displayed.
- Select the application you want to assign to the new task.
- Touch Done.
- Enter a name and a description for the task and then touch **Save**.
- The new task is saved with the factory settings and displayed in the task list.

Sorting the Task List

New tasks are always added to the end of the task list. After you have configured several tasks, it may be helpful to sort the task list.

➤ The task selection is displayed.

Sort.

(🚞)

- Touch Sort.
- The list is sorted alphabetically (A Z).
- To reverse the sorting, re-touch **Sort**.
- To sort tasks by the date of last use, re-touch Sort.

Creating New Tasks (Configuration)

The Cubis software contains the following applications that you can configure as a task to meet your own requirements:


- Weighing - Formulation
- Mass unit conversion - Weighing in percent
- SQmin Minimum sample quantity (optional)
 - Timer-controlled functions - Totalizing
- Density determination - DKD measurement uncertainty (optional)
 - Second tare (preset tare)
- Calculation - Parts counting - Checkweighing


- Individual identifiers

- Statistics
- Averaging

The basic process for task configuration is as follows:

- 1. Create new task.
- 2. Select application.
- **3.** Review all of the following prompts and configure your own settings as required.
- **4.** Perform steps 2 and 3 for additional applications if you would like to combine multiple applications.
- **5.** Check settings for weighing and printouts and modify as required.
- **6.** As Administrator, you define whether the new task can be used by all or only by certain users.
- 7. Enter a name and description for the new task.
- **8.** Save task.

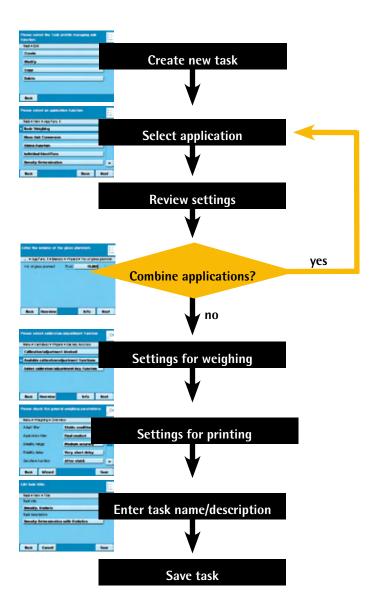

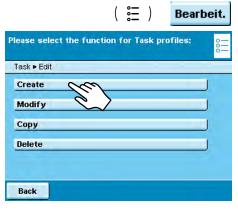


Example: Creating a task

- Open Task Management and touch Edit..
- ➤ The list of options is displayed.
- ► To define a new task, touch **Create**.



- > An application list is displayed.
- To view all applications, use the scroll bar on the right.
- ▶ Select the application you want to assign to the new task.
- ► Configure this application as required. The program guides you through the configuration menu.
- ▶ If you want to assign additional applications to the task, these applications should also be configured (see Combining Applications into One Task).

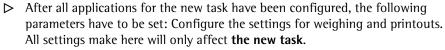




On balances verified for use as legal measuring instruments, modification of some settings may limited or prohibited altogether.

Settings for Weighing

Check all settings and change as required.

Explanations for the individual options can be found in the chapter **System Settings** in the section Device Parameters.

Touch Next.

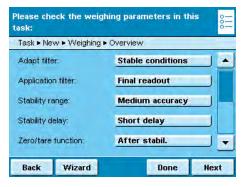

Task ➤ New ➤ Print Fct. ➤ Overview Printout interface: Com A; Com B Protocol (Com A): Print Stand, printout (Com A): GLP printing (Com A): Off Print event (Com A): PRINT

Done

Please check the printout parameters of this

Wizard

Settings for Printouts

► Check all settings and change as required.

Explanations for individual options can be found in the **System Settings** chapter in the Device Parameters section.

- Touch Next.
- Enter a short name and a description for the new task.
- Touch Save.
- ➤ The task selection list with the new task is displayed.

Configuring a Printout

You can configure the printout at two places; for example you can have the printout for basic weighing differ from all the other applications.

In System Settings (Menu)

In this menu, you define the basic settings. Previously created tasks will not be affected by modifications to the basic settings.

Creating Tasks

This is where you can specifically modify the user-defined settings for any task. You will be routinely prompted for the printer settings with every new task you create and can be modify them.

Printout (example):	Print events:	Printout content (record)	Print elements				
	Start task	21.02.2011 09:17 SARTORIUS Mod. MSA 224S SerNr. 12-34-56	GLP header				
	Start task	Laboratory technician	Operator name				
	In this example: Start task	Product	ID1, identification: fixed/variable				
	(🖹) key	123.5678g	Measured Value				
	Exit task	21.02.2010 09:20 Name:	GLP footer				



Interfaces for Printer Output

Setting Options for Printouts

The options for printouts depend on the how the interfaces and their operating modes are configured. So if not all options are available, check the interface configuration.

The configuration of standard printout parameters defines the following: Peripheral ports (Com A), US/PC (Com B), optional interface (Com C), Ethernet

(Com D)

Protocol (Operating Mode) Depends on the interface: Print, SBI, xBPI, SICS

Output Depends on the operating mode Single printout or automatic (not with printer),

with/without stability

Printout Criterion Manual without stability, manual with stability

Format Autoprint 22 characters for applications, 16 characters for raw data

Standard Data Output Standard 1 (measured value only), Standard 2 (measured value and date/time),

Standard 3 (date/time, weight block (N,T,G) and dotted line

GLP-compliant Printout Turned off/on for Cal./adjust./Always On

Print Event Defines which event should trigger a print job.

The selectable print events depend on the selected application.

For basic weighing: () key, start or end task, calibration/adjustment result;

for calculating applications: Additionally, initialization and result For logging applications: Additionally, component and evaluation


Print Elements The elements to be printed also depend on the selected application. You define the


print elements for each each event, e. g. date and time, dotted line, measured value

and the sequence they should be printed in, etc.

Tare after Printing: Turn off, on (auto. tare after printing)



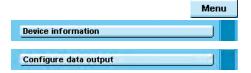

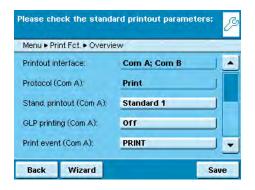

Configuring Printouts in the System Settings (Menu).


The settings can only be modified as long as no task has been activated (ready to start) in the operating mode. Make sure that basic weighing is active in the operating mode.

- Toggle to the System Settings menu.
- Open the Configure data output submenu.

If the submenu is not selectable, check the operating mode and select basic weighing.

The setting menu will open. Several options are not modifiable because they are pre-set by the interface configuration.

The initial default setting and several exemplary settings are described in the following.

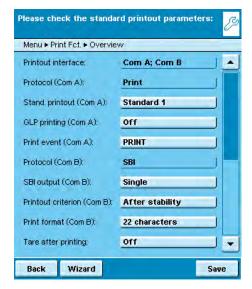

Standard printout parameters when the balance is first delivered

When the balance is first delivered, two interfaces are configured: The peripheral port (Com A) is configured for one printout, the US interface (Com B) is configured for data with a PC per SBI protocol.

Interfaces: Com A (peripherals port) and Com B (US/PC)

Protocol for Com A: Print (pre-set by the interface configuration)

Standard for output to Com A: Standard 1 (measured value only)

GLP-compliant printout to Com A: Off

Print event for output to Com A: Press the PRINT (🔁)key

Element to be output via Com A: Measured value

Protocol to Com B: SBI (pre-set by the interface configuration)

SBI output to Com B: Single printout to the PC

Criterion for output to Com B: After stability of balance

Format for output to Com B: 22 characters

Tare after printing: Off

➤ You can select the variable parameters individually by clicking on the respective input field.

or

Wizard

► To be guided step-by-step through the settings, use the wizard.

Example:


Manual printout to peripherals port, SBI operating mode

➤ After manual operation, you want an individual printout to be printed to the ported peripheral device.

Interface: Com A (peripherals port)

Protocol: SBI (toggling to the interface configuration possible)

SBI Printout: Individual printout Criterion: Manual without stability Format: 22 characters (for applications) Tare after printing: Turned on or off

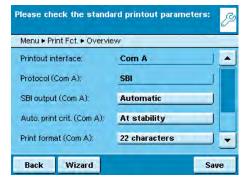

Please check the standard printout parameters:

Stop auto. print (Com A): Not possible

Auto.print interv. (Com A): 1 disp.upd.

Wizard

Automatic

At stability

22 characters

Menu ► Print Fct. ► Overview
SBI output (Com A):

Auto, print crit, (Com A):

Print format (Com A):

Example:

.

J

Save

Automatic printout to peripherals port, SBI operating mode

> At a pre-set interval, automatic printouts should be sent to the ported peripheral device.

Interface: Com A (peripherals port)

Protocol: SBI (toggling to the interface configuration possible)

SBI Printout: Automatic printout

Criterion: Without stability, at stability or after load change

Format: 22 characters (for applications)

Stop automatic printing: Not possible or by using the Print (🖹)key.

Interval for automatic printing: Number of cycles (1, 2, 5, 10 to 100) and 0.1, 0.2,

0.5 to 10 seconds.

Example:

Automatic printout to peripherals port; remote display operating mode

Please check the standard printout parameters:

Menu • Print Fct. • Overview

Printout interface:

Com A; Com B

Protocol (Com A):

Second display

Auto. print crit. (Com A):

Without stability

Print format (Com A):

22 characters

Auto.print interv. (Com A):

1 disp.upd.

At a pre-set interval, automatic printouts should be sent to the ported peripheral device as remote display.

Interface: Com A (peripherals port)

Protocol: Remote display (toggling to the interface configuration possible)

SBI Printout: Automatic printout

Criterion: Without stability, at stability or after load change


Format: 22 characters (for applications)

Interval for automatic printing: Display updates (1, 2, 5, 10 to 100) or every 0.1, 0.2,

0.5 to 10 seconds.

Example:

Printout to peripherals port only by using the PRINT (🔁) key; printer operating mode

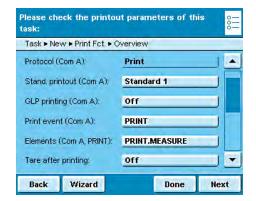

After pressing the Print key, the measured value should be printed to the ported peripheral device.

Interface: Com A (peripherals port)

Protocol: Print (toggling to the interface configuration possible)

Standard data output: Standard 1 (measured value only), Standard 2 (measured value and date/time), Standard 3 (measured value with date/time and weight block)

GLP-compliant printout Turned on or off

Print event: Press the PRINT key

Elements for the Print event PRINT: Measured value PRINT.MEASURE

Tare after printing: Turned on or off

Setting Options of Printout Parameters

Please select the standard printout (Com A): Menu > Print Fct. > Wizard > Stand. printout (Com A) > Standard 1 (measured value only) Standard 2 (measured value and date/time) Standard 3 (...) User input Back Overview Next

Standard Data Output


You can select which data should be printed out as default.

Standard 1 (measured value only)

Standard 2 (measured value with date/time)

Standard 3 (measured value with date/time and weight block)

User input is only selectable if the standard settings have been modified by the user.

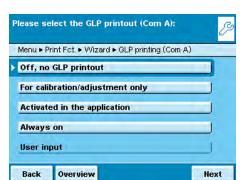

GLP-compliant Printout

You can select if and when the printout should be GLP-compliant


Off, no GLP printout

For calibration/adjustment only

Activated in the application

Alwayson

User input is only selectable as long as none of the three options apply: i.e. if only a GLP header but not a GLP footer has been configured.

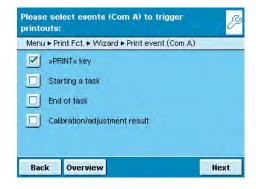

Print events

You can define one or several events that trigger a printout. The number of selectable print events depends on the selected application.

Basic Weighing:

Press the PRINT key

Starting a task

End of task

Calibration/adjustment result, after calibration/adjustment

For **Calculating applications** the following is **additionally** possible:

Initialization


Result

For **logging applications** the following is **additionally**:

Component

Evaluation

The application-specific print events cannot be configured in the Basic menu but only when you create a new task.

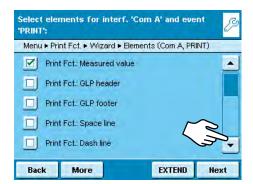

Print Elements

Which information should be printed?

Measured value

GLP header

GLP footer

Empty line

Dotted line

Form feed Date and time

etc.

- To scroll through the extensive list, slide the right scroll bar downwards.
- ► To view the complete list, touch **EXTEND**.

Reduce


Duplic

Down

Up

EXTEND

- To minimize the list, touch Reduce.
- More To
- To reset to default, touch Standard.

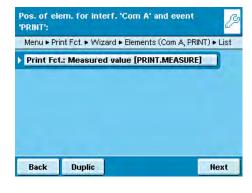


Sequence of Print Elements

In which order should the print elements be printed?

The list of selected print elements is displayed.

You can change the order and can duplicate individual elements, for multiple printouts (e. g. at the beginning and end of the list).

- Mark the element that you wish to move or duplicate.
- ► To duplicate the element, touch **Duplic**.
- ► To move the element downwards in the list, touch **Down**.
- ► To move the element upwards in the list, touch **Up**.

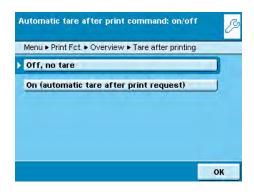

Automatic Tare after Printing

Should the balance be tared automatically after each printout?

Off, no tare

On (automatic tare after print request




Setting Options for SBI Output

SBI Output Mode

How should SBI printout be output?

Automatic print: The printout is generated automatically in the weighing cycle

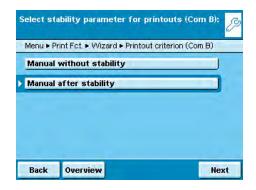


Stability Parameter for Printout

Should the printout be dependent on balance stability?

Manual without stability: Immediate printout without waiting.

Manual after stability: Printout generated immediately at stability, otherwise not until after next stability.

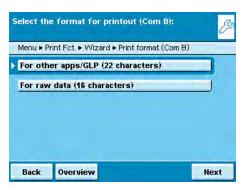

Format for Printout

How should the printout be formatted?

For other applications (22 characters): Printout with header; 20 characters plus line break (CR) and dotted line (LF)

For raw data (16 characters): Printout without header; 14 characters plus line break (CR) and dotted line (LF)

Combining Applications into one Task

You can combine applications that are saved in the balance into one task. The following rules need to be followed (see table below):

- An application can only be combined with applications from **different groups**.
- You can only combine the applications of the first three groups in the given order, e.g. weighing with checkweighing and totalizing. The sequence weighing, totalizing and checkweighing is not possible.
- The applications of the Additional Functions group can be combined freely, the order is not relevant in this case.

When you combine applications, the Cubis software automatically suggests only those applications that follow the corresponding rules.

Application group	Application	Weighing	Mass unit conversion	SOmin function	Individual identifiers	Density determination	Statistics	Calculation	Averaging	Formulation	Weighing in percent	Timer-controlled actions	Totalizing	DKD measurement uncertainty	Second tare memory	Parts counting	Checkweighing	lonizer
	Weighing	0	•	•	•	•	•	•	•	•	•	•	•	•	•	•	•	•
ng Suc	Density determination	•	•	X	•	0	•	0	0	0	0	0	0	X	•	0	0	•
Calculating applications	Calculation	•	•	X	•	0	•	0	0	0	0	•	•	X	•	0	•	•
Calc	Averaging	•	•	•	•	0	•	O	0	O	0	O	•	•	•	0	•	•
	Weighing in percent	•	•	X	•	0	•	0	0	•	0	0	•	×	•	O	•	•
	Parts counting	•	•	×	•	0	•	0	0	0	0	O		X	•	0	•	•
Check- ing applica- tions	Timer-controlled actions	•		•	•	O	•	•	O	0	•	O	0	•	•	O	0	•
Check- ing applica- tions	Checkweighing	•	•	•	•	O	•	•	•	O	•	O	•	•	•	•	O	•
	Statistics	•	•	•	•	•	0	•	•	0	•	•	0	•	•	•	•	•
Logging applications	Formulation	•	•	•	•	O	O	O	O	0	•	O	O	•	•	O	C	•
Lo	Totalizing	•	•	•	•	O	0	•	•	O	•	O	O	•	0	•	•	•
S	Mass unit conversion	•	C	•	•	•	•	•	•	•	•	•	•	•	•	•	•	•
Additional functions	SQmin function	•	•	O	•	⊼	•	⊼	⊽	•	₹	•	•	•	•	₹	•	•
fun	Individual Identifiers	•	•	•	0	•	•	•	•	•	•	•	•	•	•	•	•	•
ional	DKD measurement uncertainty	•	•	•	•	X	•	X	₹	•	⊼	•	•	O	•	×	•	•
dditi	Second tare memory	•	•	•	•	•	•	•	•	•	•	•	•	•	O	•	•	•
<	lonizer	•	•	•	•	•	•	•	•	•	•	•	•	•	•	•	•	O

- Weighing is always available
- Combination possible
- Combination possible is only activated as long as no calculating application is active.
- No combination with identical application possible
- No combination with an application of the same group possible
- No combination possible because it is not meaningful

Executing Tasks

After all required settings have been configured, you can now start and execute a task.



- ▶ If you haven't already done so, go to Task Management.
 - > The task selection is displayed.
- Touch the desired task.





- ▶ If the desired task is already selected (dark background and arrow on the left), touch Start.
- > The program switches to operating mode and the selected task is activated.
- ► Follow the instructions on the display.

Weighing


Purpose: Weight determination within the scope of the device-specific weighing capacity

(see Specifications).

Combination options: Mass unit conversion, second tare memory and individual identifiers,

(SQmin function, DKD measurement uncertainty).

Configure Task: My Weighing

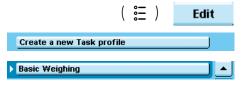

Stable conditions

Medium accuracy

Final readout

Short delay

Please check the weighing parameters in this

Task ► Weigh ► Weighing ► Overview

Adapt filter:

Application filter:

Stability range:

Stability delay:

Back

Cancel

Weighing without combining it another application:

- The settings for Weighing are displayed.
- Check all settings and change as required.
- Touch Next.
- ▶ The settings for **Printouts** are displayed.
- ▶ Check all settings and change as required.
- Touch Next.
- ▶ If you are administrator: You define whether all or only individual users may use this task.
- Touch Next.
- ▶ Enter a short name and a description for the new task (e. g. "My Weighing").
- To save the new task, touch Saue.



Execute Task: My Weighing

- ▶ If you haven't already done so, go to Application Management.
- ➤ The task selection is displayed.
- ► Touch the desired task.

Start

or

Save

- ▶ If the desired task is already selected, press **Start**.
- ▶ The program switches to the operating mode, and the task is activated.





Opening and closing the rotation draft shield (models MSx6.6S/3.6P/2.7S):




Defining the opening width:

Touch DraftShd.

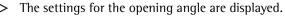

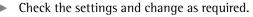

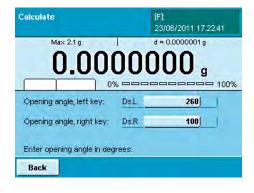



Draft shield opens according to the opening angle setting.

Mass Unit Conversion

Purpose: The weight value can be displayed in up to a maximum of five different units in

succession (see table below). However, the desired weight units must be assigned

beforehand (1 through 5).

Combination options: Second tare, (SQmin function, DKD measurement uncertainty)




Not all units programs can be selected on balances used as legal measuring instruments. The user can only switch between metric units (e. g. milligram, gram, kilogram) (also model-specific).

Conversion Factors for Weight Units

Unit	Factor	Display				
Gram	1,0000000000	g				
Kilogram	0,0010000000	kg				
Carats	5,0000000000	ct				
Pounds	0,00220462260	lb				
Ounce	0,03527396200	OZ				
Troy ounce	0,03215074700	ozt				
Hong Kong taels	0,02671725000	tlh				
Singapore taels	0,02645544638	tls				
Taiwan taels	0,0266666000	tlt				
Grains	15,43235835000	GN				
Pennyweight	0,64301493100	dwt				
Milligram	1000,0000000000	mg				
Parts per pound	1,12876677120	/lb				
Chinese taels	0,02645547175	tlc				
Mommes	0,26670000000	mom				
Carats	5,0000000000	ct				
Tola	0,08573333810	tol				
Baht	0,06578947437	bat				
Mesghal	0,21700000000	MS				
Ton	0,00000100000	t				
Pound:Ounce	0,03527396200	lb:oz				
Newton	0,00980665000	N				
Microgram	1000000,00000000000	μg				


Configure Task: Mass Unit Conversion

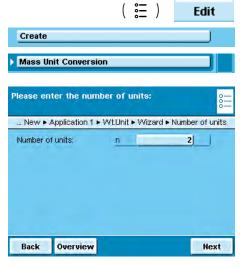

► Select the desired number of units (max. 5) to be used to display the weighing results.



- Now you will be prompted to enter the display accuracy and weight units of all the desired units one by one.
- Select the **Weight unit** (e. g. gram, kilogram, carat, pound) for the first unit.
- Select the **Display accuracy**: (e. g. complete display of all decimal places or leave off last place).

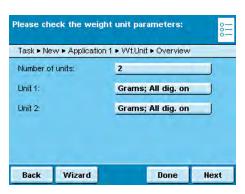

- ➤ An overview of all settings is displayed.
- Check settings and change as required.
- ➤ A message asks whether you would like to add another application.
- ▶ If required, select additional applications to be combined.
- Now you will be prompted to configure the weighing and printing functions.
- Check all settings and change as required.
- Enter a short name and a description for the new task.
- ► To save the new task, touch **Save**.

Execute Task: Mass Unit Conversion

(등)

- ▶ If you haven't already done so, go to Application Management.
- ➤ The task selection is displayed.
- ► Touch the desired task.

or

Start

- If the desired task is already selected, press **Start**.
- Follow the instructions on the display.



Unit

- ► To change the displayed unit, touch **Unit**.
- ➤ Select the unit to be used to display the weighing results.

 This setting remains in effect until you change the selection or end the task.
- ► To display weighing results in all units, touch **Toggle**.



Minimum Sample Quantity SQmin

Purpose:

This application is used to compare the measured weight value directly with the defined minimum sample quantity (SQmin = sample quantity minimum). This ensures that weighing results are within tolerances defined by your quality assurance system. This fulfils the requirements of United States Pharmacopeia (USP). According to USP guidelines, the measurement uncertainly can be a max. of 0.1% related to the initial weight when weighing substances for quantitative analyses.

Requirements:


The balance must be set up by a service technician to be able to use the SQmin function. The technician will determine the permitted minimum sample quantity and load this to your balance using the guidelines of your QA system. He will document this setting via a "Weighing module test as per USP" certificate in which the measurements and min. sample quantity are logged. The SQmin function ensures that the weighing results correspond to USP guidelines.

Combination options:

Mass unit conversion, second tare memory, individual identifiers, (DKD measurement uncertainty)

Edit

Configuring Task: Minimum Sample Quantity SQmin

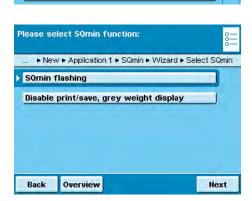

Create

SQmin Function

► Select which function should be triggered for minimum sample quantity. **SQmin flashing**: The SOmin tab flashes.

Disable print/save...: Printout and saving of the weight value are disabled; the weight value is displayed in gray.

- > An overview of all settings is displayed.
- Check all settings and change as required.
- Combine the task profile with another application?

 Task New Application 2

 No further application function

 Mass Unit Conversion
 Individual Identifiers

 DKD Uncert. of Measuremt.

 Second Tare (Preset Tare)

 Back

 Done Next
- > A message asks whether you would like to add another application.
- ▶ If required, select additional applications you would like to combine.
- > Now you will be prompted to configure the weighing and printing functions.
- ► Check all settings and change as required.
- ▶ Enter a short name and a description for the new task.
- To save the new task, touch Save.

Execute Task: Minimum Sample Quantity SQmin



- ▶ If you haven't already done so, go to Application Management.
- ➤ The task selection is displayed.
- ► Touch the desired task.

or

Start

If the desired task is already selected, press **Start**.



- ▶ When you start a task with the SQmin function, the defined SQmin value is shown on the display.
- ➤ As long as the weight value is lower than the SQmin value, the tab "SQmin" will flash.

Individual Identifiers

Purpose:

You can define identifiers for the following print jobs:

- Printout by pressing PRINT
- Task start
- Task end
- Initialization (for calculating applications only)
- Result (for calculating applications only)
- Component (for logging applications only)
- Evaluation (for logging applications only)

These identifiers are also printed in the log.


Combination options:


Second tare, (SQmin function, DKD measurement uncertainty)

Example:

You want to configure the application as follows:

- You want the printout to be activated with the PRINT key.
- You want to configure the application so that four identifiers are included in the printed log, i.e. your company address in three lines and the batch number in one line
 - The batch number is to be counted up automatically.

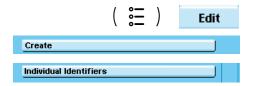


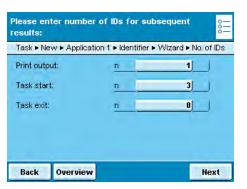


▶ Enter the desired number of individual IDs that you want to set up.

Print output: Identifier for the Print key

Task start, task end: Number of identifiers for task start or task end.

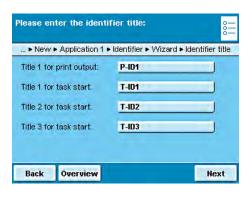

- > The corresponding ID (P-ID1, T-ID1, T-ID2, etc.) will be displayed for each identifier.
- Enter the desired title for each ID.

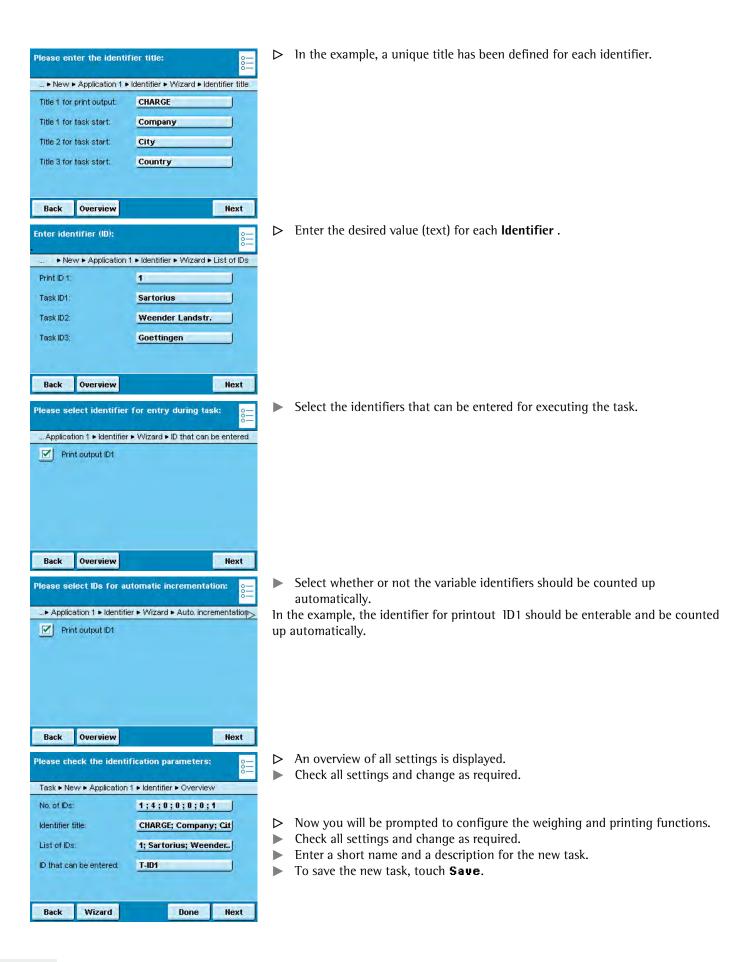


Execute Task: Individual Identifiers

- ▶ If you haven't already done so, go to Application Management.
- ➤ The task selection is displayed.
- ► Touch the desired task.

or

Start

▶ If the desired task is already selected, press **Start**.

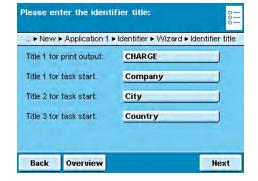

> The defined identifiers are displayed (cf. example above).

Touch Next.

Now the predefined lines are printed when the task starts:

Firma: Sartorius Strasse: Weender Landstr. Ort: Goettingen



- Place the sample on the weighing pan and press the PRINT key \Box .
- ➤ The identifier for the printout is displayed and can be changed, if necessary.
- ▶ In the example, CHARGE was defined for the printout.
- ▶ If required, you can change this value now.
- ► Touch **Next** or the PRINT key 🖯 once more.



- ➤ To log the value, press the Print key.
- ➤ The complete printout looks like this:

Firma: Sartorius
Strasse:
Weender Landstr.
Ort: Goettingen
N + 214.5 g
CHARGE 1
N + 213.3 g
CHARGE 2

Density Determination

Purpose: You can determine the density and volume of liquid, solid or pasty materials.

Combination options: Checkweighing, statistics, second tare or individual identifiers.

Requirements: You require the Sartorius YDK01MS density determination kit or your own

customer-specific set. The basic calculations for density determination are described

at the end of this chapter.

Features *= Factory setting

Liquid

*Water: for density determination using water; the density for the corresponding

temperature is calculated automatically

Ethanol: for density determination using ethanol; the density for the corresponding

temperature is calculated automatically

Individual entry: for density determination using another liquid, enter liquid and its density for

corresponding temperature

Density at 20°C and coefficient: for density determination using another liquid, enter density at 20 °C and the

coefficients of expansion, the density for the corresponding temperature

is calculated automatically.

Air density

*Air density at 20°C: for measurement under standard laboratory conditions

Individual entry: for measurement under other conditions, enter air density

Decimal places

None: The measurement value is displayed without decimal places.

1, *2, 3 or 4 places: The measurement value is displayed with the selected number of decimal places.


You can choose from **four methods** for density determination:

- Determine density of liquid (with glass plummet)


- Determine density via buoyancy with Sartorius YDK01MS density determination kit (for solid samples)
- Determine density via displacement (for solid samples)
- Determine density via pycnometer (for liquid, pasty and powder samples)

Configure Task: Determining the Density of a Solid

You can choose from two methods to determine the density of a solid: **Buoyancy** or displacement.

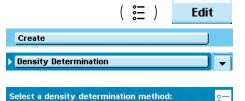

Select the method that you want to use: Buoyancy or displacement.

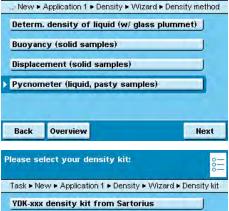

- Select the density kit you want to use.
 - Sartorius density kit YDK: For measuring with the YDK kit A different density determination kit: If you want to use another kit No density kit: If you do not want to use any kit.
- Select or enter the prompted values (for explanations see **Features** on page 78).



Additional prompts/settings for another determination kit


*= Factory setting

Parameters for the density kit

Name: Enter a name


No. of wires: Enter the number of wires for the density kit that are immersed into the liquid. Wire diameter: Enter wire diameter (in mm)


Vessel diameter: Enter vessel diameter (in mm)

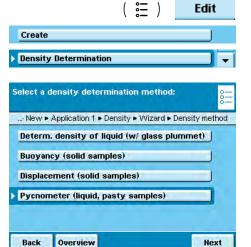

Configure Task: Determining Density Using a Pycnometer

- Select the **Pycnometer** method.
- Select or enter the prompted values (for explanations see Features on page 78).

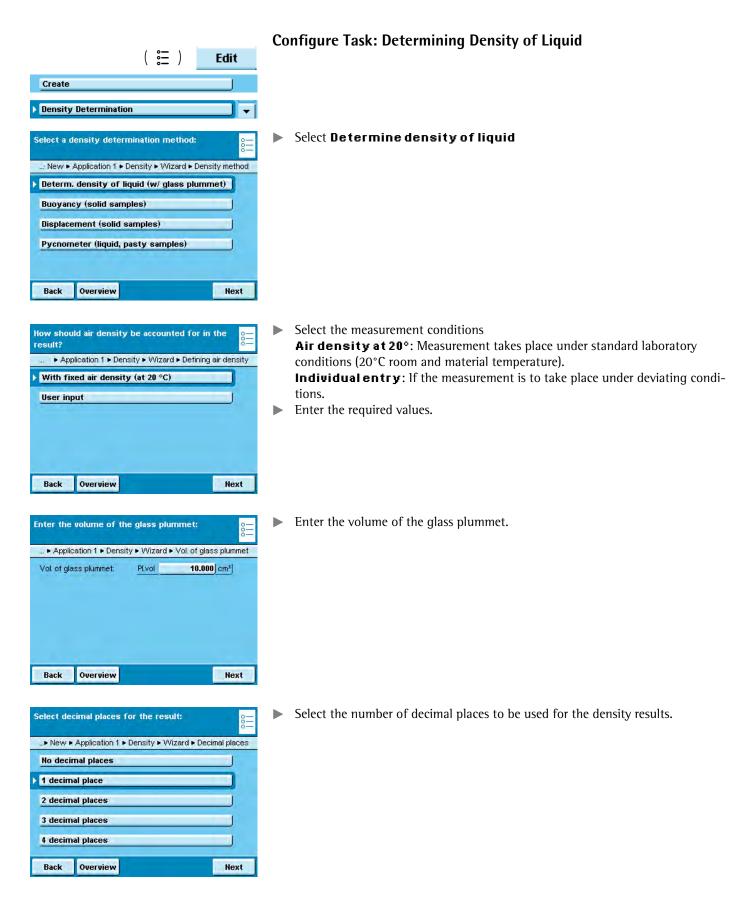

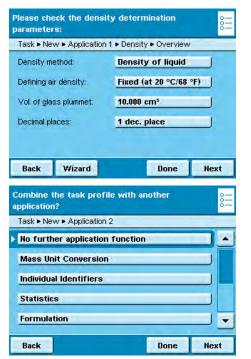

Check all settings in the overview and change individual parameters as required.

- ➤ An overview of all settings is displayed.
- Check all settings and change as required.
- ➤ A message asks whether you would like to add another application.
- ▶ If required, select additional applications to be combined.
- Now you will be prompted to configure the weighing and printing functions.
- Check all settings and change as required.
- Enter a short name and a description for the new task.
- To save the new task, touch Save.

Execute Task: Density Determination



- Prepare the density determination kit that you want to use
- ▶ If you haven't already done so, go to Application Management.
- > The task selection is displayed.
- ► Touch the desired task.

or

Start

▶ If the desired task is already selected, press **Start**.



- > The program switches to operating mode.
- ► Follow the instructions on the display.

Calculation Basis for Density Determination

The density determination application is based on the equations:

RhoDensity of sample (ρ) $ρ_{fl}$ Density of buoyancy liquid

ρ_{fl} Density of buoyancy liquW_a Weight of sample in air

Wfl Weight of sample in liquid (for determining density of liquids, buoyancy and displacement)

Weight of fill medium (for pycnometer)

W_r Weight of sample and fill medium (for pycnometer)

LA Air buoyancy correction = 0.0012 g/ccm

Buoyancy: Rho = $(W_a * (\rho_{fl} - LA)) / ((W_a - W_{fl}) * corr) + LA$

Displacement: Rho = $(W_a * (\rho_{fl} - LA)) / (W_{fl} * corr) + LA$

When the displacement method is used to correct for buoyancy by means of a wire (string) suspended in liquid, the factor 1.00000 is used in the calculation as factory-set default.

The correction factor **Corr** is calculated from: $Corr = 1 - n * d^2 / D^2$ with:

n No. of wires:

d Wire diameter (YKD01MS: 0,7 mm)

D Inner diameter of the vessel being used (YKD01MS: 76 mm)

Please check the density kit parameters:

Task ▶ New ▶ App.Func. 1 ▶ Density ▶ Wizard ▶ D.kit param.

Name of kit:

Density set 2

Number of wires:

n 2

Wire diameter:

d 1.00 mm

Vessel diameter:

D 76 mm

When other vessels or other submersion assemblies are used, this factor can be changed by modifying the parameters for the density kit.

Statistics

Purpose:

Saving and statistical analysis of weight values and other calculated values. The values are generated as results:

- Transaction counter
- Sum of all values
- Average value
 - Standard deviation
 - Lowest value (minimum)
 - Highest value (maximum)
 - Difference between maximum and minimum

Options:

You can save a maximum of 65535 components.

Combination options:

Density determination, calculation, averaging, weighing in percent, parts counting, timer-controlled functions, checkweighing, mass unit conversion, second tare memory, individual identifiers (SQmin function, DKD measurement uncertainty)

Create Edit

Configure Task: Statistics


- * = factory setting
- ▶ Define how the weight value will be applied.
- *Manually: The weight value is applied as soon as the corresponding button is touched.

Automatically: first value at stability: The weight value is applied automatically as soon as the balance has stabilized after filling and the value exceeds the minimum load.

Automatically: last value at stability: The last weight value with stability is applied automatically, as soon as the balance is unloaded and the value falls short of the minimum load. If stability was reached repeatedly between measurements, the last stable weight value will be stored.

You must now select the minimum load to be applied:

(none, 10, 20, ...*100 ... 1000 digits).

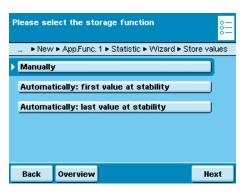



Weighing Mode with Automatic Storage:

▶ Select the weighing mode to be used for the weight values.

Initial weighing: Initial weighing of samples in a container.

Final readout: Final weighing of samples in a container.

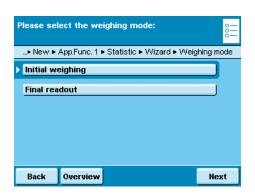

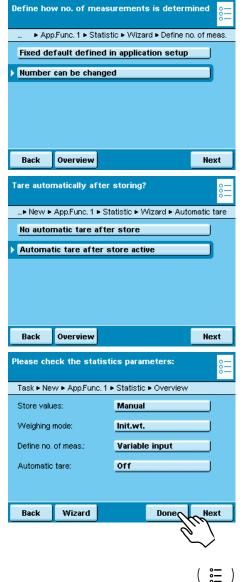

Define when the number of measurements (items) will be defined.

Enter fixed number: The item number is defined beforehand.

Number can be changed: Enter the item number when you execute the task.

▶ Define whether or not automatic taring takes place after the transfer.

- An overview of all settings is displayed.
- Check all settings and change as required.
- ▶ A message asks whether you would like to add another application.
- ▶ If required, select additional applications to be combined.
- Now you will be prompted to configure the weighing and printing functions.
- Check all settings and change as required.
- Enter a short name and a description for the new task.
- To save the new task, touch Save.
- ▶ The new task is saved with your settings and displayed in the task list.

Execute Task: Statistics

- If you haven't already done so, go to Task Management.
- ➤ The task selection is displayed.
- ► Touch the desired task.


or

Start

If the desired task is already selected, press Start.


▶ The program switches to operating mode

- ▶ If required, enter the number of items that you want to analyze.
- ▶ Place the first sample on the weighing pan.
- ► To start Statistics, touch Start.
- > The first weight value is displayed.

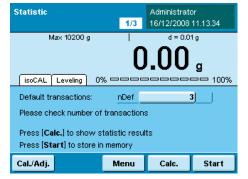



- ▶ Place the next sample on the weighing pan and touch **Next**.
- ► Continue with all other samples in the same way.
- You can remove each last value from the statistics by touching Correction.
- During the task, you can have the result displayed at any time by touching **Calc.**.

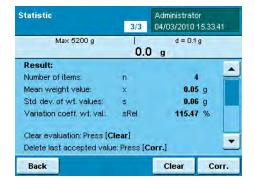

- > Once all intended samples have been weighed, the result is displayed.
- ► To delete the results and begin the next task, touch **Clear**.

Print example:

n	1	
Comp	98.8	g
n	2	
Comp	56.4	g
:		
:		
:		
n	5	
Comp	56.4	g
n	5	
Sum	280.3	g
X	56.06	g
S	24.93	g
sRel	44.46	%
Max	98.8	g
Min	37.2	g
Diff	61.6	g

Calculation

Purpose: This application is used to calculate the weight value using an algebraic formula

that you define yourself. You can define a fixed formula or one that you can


change while executing the task.

On balances used as legal measuring instruments, follow the directions in the "Executing tasks" section!

Combination options: Second tare, (SQmin function, DKD measurement uncertainty)

Example: You want to determine the area of irregularly cut pieces of paper. The grammage

of the respective paper type is known.

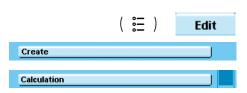

...New ▶ App.Func. 1 ▶ Calculate ▶ Wizard ▶ Defining formula


Define how the formula is determined

Fixed formula

Back

Variable formula (input)

Overview

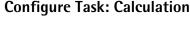

➤ Select whether you want to use a fixed formula or a variable formula.

Enter Fixed formula: All calculations are carried out with the formula that you

Formula can be changed: You can change the formula before each calculation.

Example: The area of irregular paper samples must be determined.

The respective paper grammage in g/cm² is known. The equation for determining the paper area in cm² is:

Paper area = weight value W in $g * 10000 / grammage in <math>g/cm^2$.

The formula to be entered for a grammage of 80 g/cm² is:

Wx10000/80

Next


define for this task.


To analyze different paper types, the formula must be able to be changed before each calculation (e.g. for paper with 80, 90, 100 g/cm^2).

Additional formula examples are at the end of this section. Likewise, you can create several tasks, one for each paper type.



▶ If you want to work with a variable formula, enter the formula when you execute the task.

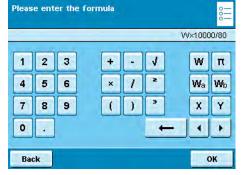


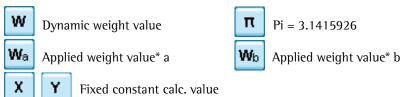


- Square root**

 2 2. Power **


 3 3. Magnitude**
- * for calculation
- ** corresponding to general mathematical characters

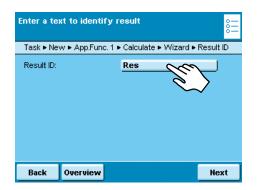


► Enter a name for the event print e. g. "Area" (max. 6 characters).

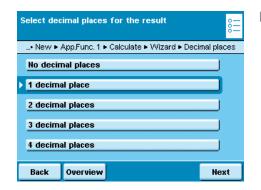

Define how many decimal points should be used for the results: none,*1...7 digits

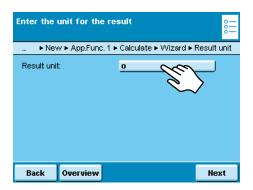

► Enter the unit to be used for the results (max. six characters, e.g. cm²)

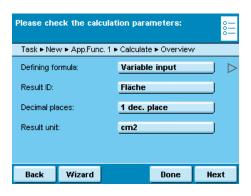

- ➤ An overview of all settings is displayed.
- Check all settings and change as required.

A message asks whether you would like to add another application.

- > Prompts for configuring weighing and printing appear.
- Check all settings and change as required.
- ► Enter a short name and a description for the new task.
- To save the new task, touch **Save**.

Examples of Calculation Formulas

1) Diameter determination for cylindrical bodies

Definition:

Diameter = $2 \times V (W / (\pi \times I \times Rho))$


with: V = Square root

W = weight value in g

 $\pi = 3.145$

l = length of sample in cm

Rho = density of sample in g/cm^3 (e.g. 8.30000 g/cm^3)

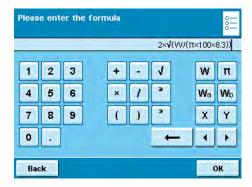

Formula

Input:: $2 \times V (W/(\pi \times 100 \times 8.3/1,000))$

Header: d Unit: mm

2) Air buoyancy correction

Purpose

This application can correct weighing errors that can arise due to air buoyancy when working with weights of different densities.

The air density value is required when calculating the air buoyancy correction.

Formulas and abbreviations

1. Definition: m = W * K

with: m = mass of the sample, unit symbol "g!"

W = weight value in grams K = correction factor


2. Definition: K = 1 - (RhoL / RhoST) / 1 - (RhoL / RhoG)

with: RhoL = air density in g/cm³, e.g. 0.0012 g/cm³ (standard air density)

RhoST = density of steel, e.g. 8.0000 g/cm³ RhoG = density of sample, e.g. 2.4000 g/cm³

Preparation

Access the "Change" item from the Task menu and select the "TASK-AIR-BUOYANCY_CORR:" task downloaded to the SD card.

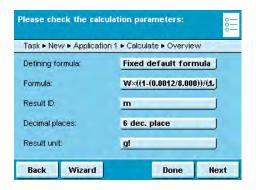

The "Calculation" application is set with the following parameters:

- Formula specifications: fixed

Formula: $W \times ((1 - (0.0012 / 8.000)) / (1 - (0.0012 / X)))$


- Result identifier: m

- Number of


decimal places: 6 digits
- Result unit: g!

Executing a Task: Calculation "Air buoyancy correction"

- ▶ If you haven't already done so, go to Task Management.
- ➤ The task selection is displayed.
- ► Touch the "Air buoyancy correction" task from the Task menu.
- ▶ The formula for the air buoyancy correction is already defined.
- ► Enter the density of the sample via "Variable X," in this example 2.400 g/cm³.
- Touch Start to begin the calculation.

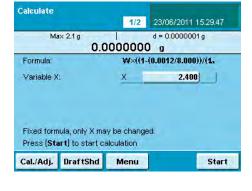

- ▶ The calculation results are displayed.
- To return to the weight value display, touch **Weigh**.
- ► You can view both values, the weight and the mass.

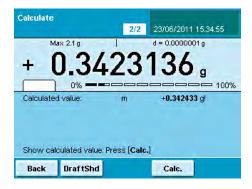

- ▶ If you want to carry out additional weighing and calculations using the same formula, place the sample on the balance and then touch Calc.
- ► If you want to carry out additional weighing using a different formula, touch **Back**.
- ▶ You can now change the formula and then proceed as described above.

```
Print via the ( \bigcirc ) key: N + 3.183629 g X 2.400 m + 3.184744 g!
```

3) Diameter determination

Purpose

This application is used to determine the diameter of round wires and metal threads (cylindrical solid bodies). For example, for use in determining the diameter of filaments. The density and length of the sample is required for calculating the diameter.

Formulas and abbreviations

Definition: $d = 2 \times \sqrt{(W / (\pi \times I \times Rho / 1000))}$

with: d = diameter of sample in mm $\sqrt{ = square root}$

W = weight value in grams

vv = weight value in grains

 $\pi = Pl = 3,145..$

l = length of sample in mm, e.g. 100 mm Rho = density of sample in g/cm³, e.g. 8,3 g/cm³

Preparation (only required if you need to make changes)


Access the "Change" item from the Task menu and select the "TASK-DIAMETER-DETERM" task downloaded to the SD card.

The "Calculation" application is set with the following parameters:

Formula specifications: fixed


- Formula: $2 \times \sqrt{W/(\pi \times X \times 8.3/1000)}$

Result identifier: dNumber of decimal places: 3 digitsResult unit: mm

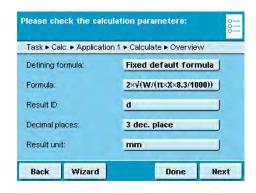

Executing a Task: Calculation "Diameter determination"

- ▶ If you haven't already done so, go to Task Management.
- ➤ The task selection is displayed.
- ► Touch the "Diameter determination" task from the Task menu.
- ➤ The formula for diameter determination and the density of the sample (8.3 g/cm³) are preset..
- ► The length of the sample, e.g. 100 mm can be entered using the X variable.
- ► Touch **Start** to begin the calculation.

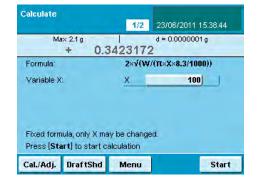

- ► The calculation results are displayed.
- To return to the weight value display, touch **Weigh**.
- You can view both values, the weight and the diameter.



- ► If you want to carry out additional weighing and calculations using the same formula, place the sample on the balance and then touch Calc..
- If you want to carry out additional weighing using a different formula, touch **Back** Back.
- ▶ You can now change the formula and then proceed as described above.


Execute Task: Calculation




- If you haven't already done so, go to Application Management.
- The task selection is displayed.
- Touch the desired task.

Start

If the desired task is already selected, touch Start.

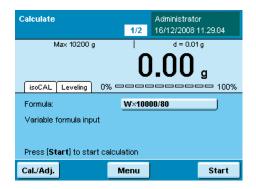

- When you want to work with a variable formula, you can now enter or change the formula.
- Place the sample on the balance.

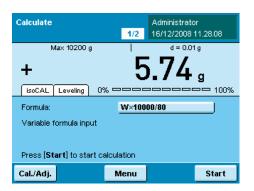


- The weight value is displayed.
- To start the calculation, touch **Start**.



- The calculation results are displayed.
- To return to the weight value display, touch Weigh.

If you want to carry out additional weighing and calculations using the same formula, place the sample on the balance and then touch Calculate. Back

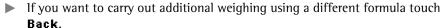

You can now change the formula and then proceed as described above.

Averaging

Purpose:

This application is used for moving samples (e. g. live animals) and for weighing in unstable environments. A measurement cycle is automatically carried out with a defined number of measurements for each object to be weighed. The individual measurements are averaged and this average is displayed as the result.



On balances used as legal measuring instruments, follow the directions in the "Executing tasks" section!

Options:

You can start the measurement manually or it can start automatically as soon as the min. weight is reached. Depending on how strong the unsteadiness of the sample is, you can vary the sensitivity for the measurement. The number of measurements from which the average is taken can also be defined.

Combination options:

Checkweighing, timer controlled functions, totalizing, statistics, formulation, SQmin function, DKD measurement uncertainty, second tare, individual identifiers.



Configure Task: Averaging

Select how the application should start.

Manual start by pressing key: This starts the application manually. *Automatic start: The application starts as soon as a pre-defined min. load is exceeded.

Select minimum load that triggers start

Task ► New ► App.Func. 1 ► Averaging ► Wizard ► Min.load

10 digits

20 digits

50 digits

► 100 digits

Back Overview Next

Please select the weighing mode:

■...New ► App.Func. 1 ► Averaging ► Wizard ► Weighing mode

Initial weighing

Final readout

Next

Back

Overview


▶ Define the min. load for the automatic start of the averaging and/or for the release of the result display for the manual start.

none, 10, 20, 50, *100, ... 1000 digits

The number of **digits** depends on the balance resolution (see Display). Example: For a balance with a resolution of d = 0.1 (100 mg), 100 digits means a min. load of $100 \times 0.1 = 10$ g. With this setting, the automatic start would be triggered as soon as a min. of 10 g was placed on the balance.

Only for automatic start:

Determine the weighing mode: *Initial weighing or Final readout.

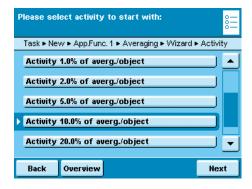


► Enter the range of the unrest for the start of the measurement: 0.1%, 0.2%, 0.5% ... *10% ... 50% from measured object


This entry estimates the strength of unrest that is caused by the acti

This entry estimates the strength of unrest that is caused by the activity (e. g. animal movement). You can use the rough estimation or a more detailed estimation using percentages.

The weight value changes with each movement of the weighing pan. As soon as three consecutive measurement values are within the selected unrest range, the interval measurement begins.

Example: You are weighing animals with an average weight of approx. 300 g. With the "5% from measured object" setting, the measurement starts when the movements are within a range of 15 g, i.e. \pm 7.5 g. With the "50% from measured object" setting, the interval measurement would start when the weight value was within a range of 150 g, i.e. \pm 75 g.

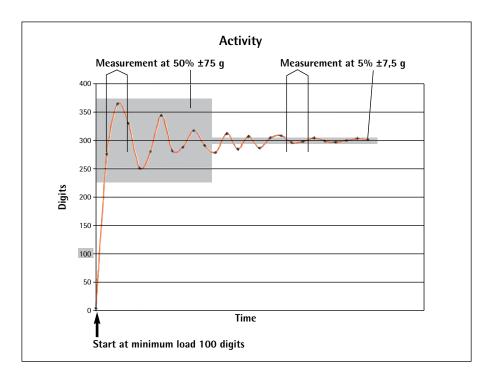


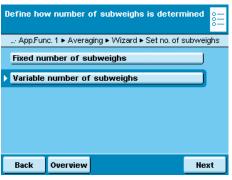

▶ Define the number of measurements to be used for the average value.

Enter fixed number: You must now enter a fixed number.

*Number can be changed: You will define the number when you execute the task.

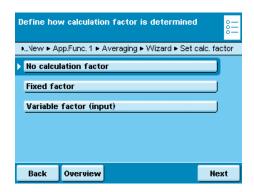


▶ Define whether or not the weighing results should be calculated with an additional factor and how this factor should be applied.

No calculation factor

Fixed factor: You must now enter this factor.

Variable factor. You must enter this factor when you execute the task.


Example: Each animal should receive a food supplement of 50 mg per kg of body weight. For this, the factor 0.05 can be entered (weight value in $g/1000 \text{ g} \times 0.05 \text{ g}$). During the measurement, both the weight value (animal weight, e.g. 285 g) as well as the calculated value (amount of supplement, e.g. can be displayed.

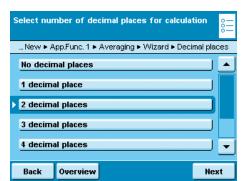

If you selected the calculation:


▶ Define the number of decimal places to be used for the calculation.

none, 1, *2, ... 6 digits

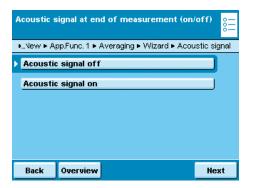

▶ Define whether or not the end of the measurement (and if necessary the calculation) should have an acoustical signal.

*Acoustic signal off or Acoustic signal on

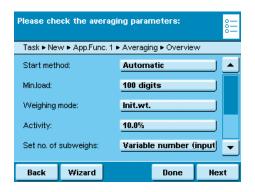

- > An overview of all settings is displayed.
- Check all settings and change as required.
- ➤ A message asks whether you would like to add another application.
- ▶ If required, select additional applications to be combined.
- > Prompts for configuring weighing and printing appear.
- Check all settings and change as required.
- Enter a short name and a description for the new task.
- To save the new task, touch **Save**.

Execute Task: Averaging with Manual Start


(📜)

- ▶ If you haven't already done so, go to Application Management.
- ➤ The task selection is displayed.
- ► Touch the desired task.

or

Start


▶ If the desired task is already selected, press **Start**.

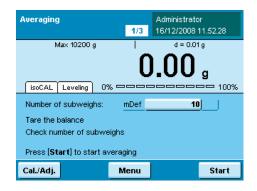

- ▶ When you are using a variable number of measurements, enter the desired number.
- ▶ When you are weighing live animals, place the container (cage) on the weighing pan and tare the balance.
- Place the animal in the container.

Start

To begin weighing, touch **Start**.

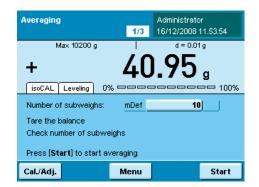


The defined number of measurements is carried out and the average is displayed.

Start

Calc.

► To begin the calculation, touch **Start**.

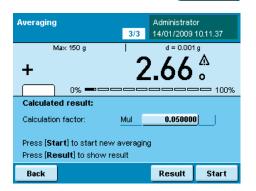

> The results calculated with the entered factor are displayed.

Result To display the average weight value, touch **Results**.

► To return to the display of the calculated value, touch Calculate..


Execute Task: Averaging with Automatic Start

- ▶ If you haven't already done so, go to Application Management.
- ➤ The task selection is displayed.
- ► Touch the desired task.

or

Start


If the desired task is already selected, press **Start**.

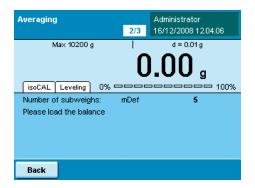


- ▶ When you are using a variable number of measurements, enter the desired number.
- ► When you are weighing live animals, place the container (cage) on the weighing pan and tare the balance.

Start

Press Start to begin the task.

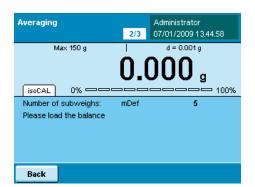

Place the animal in the container.

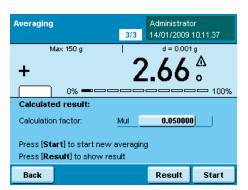

- The measurement starts automatically, the defined number of measurements and averaging is carried out.
- ➤ The calculated results are displayed.

Result

To display the average weight value, touch Results.

Calc.

To return to the display of the calculated value, touch Calculate..

Formulation

Purpose:

Combination options:

This application is used to weigh while adding several components consecutively. The balance is tared automatically after each component. The weight values of all individual components as well as the total weight results are recorded and can be logged. The weight values and the component counter are saved in protected memory so that an interrupted formulation process can be continued after the balance has been turned off or after a power failure. You can save a maximum of 65535 components.

Mass unit conversion, second tare memory and individual identifiers, (SQmin function, DKD measurement uncertainty)

Options: The program allows 7 different formulations.

1) Fixed formulation in grams: Number of components and the weight portion of each component (in g) are fixed.

2) Fixed formulation in % with total = 100%, Total weight is entered: Number of components and the percentage of each component (in g) are fixed. The total weight is entered when the task is started. If the percentages do not add up to 100, the individual values are converted to 100%.

Example: Number of components nDef =3; component 1 = 25%, component 2 = 30%, component 3 = 10%; components, total = 65% is set equal to 100%. Component 1: 25%/65% * 100 = 38.46%; Component 2: 30%/65% * 100 = 46.15%; Component 3: 10%/65% * 100 = 15,38%

- **3)** Fixed formulation in % with total = 100%, total weight is calculated according to the first component: Number of components and the percentage of each component (in g) are fixed. If the percentages do not add up to 100, the individual values are converted to 100% (cf. Formulation 2). After weighing the first component, the total weight (100%) is calculated.
- **4)** Fixed formulation in % with total <> 100%, total weight is entered: Number of components and the percentage of each component (in g) are fixed. If the total of percentages does not equal 100, the individual values are **not** converted. The total weight is entered, for each component, the user-defined weight and the percentage are displayed before weighing.
- **5) Fixed formulation in** % with total = 100%, total weight is calculated according to the first component: Number of components and the percentage of each component (in g) are fixed. If the percentages do not add up to 100, the individual values are converted to 100%. For the first component, the user-defined weight is displayed before weighing in percent. After weighing the first component, the total weight (100%) is calculated. Before weighing any other component, the respective user-defined weight is displayed in g and %.
- **6) Variable formulation in grams, component count are in a fixed order**: Before weighing the first component, you can change the number of components; before weighing each component you can change its name.
- **7) Variable formulation in grams, component count are in a fixed order**: Before weighing the first component, you can zero the number of components; before weighing each component you can change its name.

For all formulations, the following applies:

- The individual components can be stored manually or automatically (first reproducible value).
- If automatic storage of values was selected, the minimum load is determined when the task is created.
- After weighing the first component, the **Calc.** button is displayed. That way, you can change the display to show the intermediate result.
- Once the intermediate result is displayed, you can delete the value of the last weighed component by pressing the button **Corr.**; and it will be removed from the calculation. This allows you to delete all components step-by-step in reverse order. Deleting is **not undoable!**
- Use the **Back** button to toggle the display and continue the task.
- After the defined number of components has been weighed, the evaluation is displayed. Here too, you can still delete the components (beginning with the last component) and weigh them again.

Note: For the seventh formulation, there is no automatic evaluation because the defined number of components is zero. You have to operate the evaluation manually by touching Calc..

(∷) Edit Create Formulation Please select the storage function: ▶ New ▶ Application 1 ▶ Formul. ▶ Wizard ▶ Store values Manually Automatically: first value at stability Back Overview Next

▶ New ▶ Application 1 ▶ Formul. ▶ Wizard ▶ Default items

Next

Quantity can be changed during the process

Define when the no. of components is

determined

Back

Enter fixed quantity

Overview

Calc.

Corr.

Back

Configure Task: Formulation


- * = factory setting
- Define how the weight values will be applied.

*Manually: The weight value is applied as soon as the corresponding button is touched.

Automatically: first value at stability: The weight value is applied automatically as soon as the balance has stabilized. You must now select the minimum load to be applied (10, 20, ... *100 ... 1000 digits).

Set when the number of components will be defined.


Enter fixed number: The number of components is defined beforehand. *Number can be changed: Enter the number of components only when you execute the task. (Formulations 6 and 7).

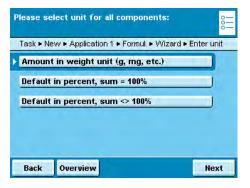

- Once you fix the components, now define now the unit:
 - *Amount in weight unit (Formulation 1)

 Default in percentage, sum = 100% (Formulations 2 and 3)

 Default in percentage, sum <> 100% (Formulations 4 and 5)
- Touch Next.
- Enter the number of components.

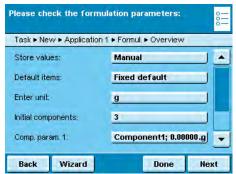

- ▶ An overview of all settings is displayed (here, Formulation 2 as an example).
- Check all settings and change as required.
- A message asks whether you would like to add another application.
- ▶ If required, select additional applications to be combined.
- Prompts for configuring weighing and printing appear.
- Check all settings and change as required.
- ▶ Enter a short name and a description for the new task.
- To save the new task, touch Save.

Execute Task: Formulation

(등)

Start


- ▶ If you haven't already done so, go to Application Management.
- ➤ The task selection is displayed.
- ► Touch the desired task.

or

▶ If the desired task is already selected, press **Start**.



- If required, enter the number of components that you want to weigh while adding them.
- To start the formulation, touch **Start**.

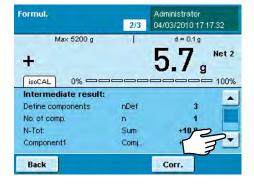

▶ If required, enter the name of components.

Please note that you will be able to scroll through the lines as soon as a scroll bar appears at the right of the display.



▶ The evaluation is displayed as soon as all components of the formulation have been added and weighed.

You can also see the complete list here when you touch the scroll bar.

Corr.

► To delete the last component, touch **Correction**.

Param.

To go to the parameter page, touch Param.

Start

To restart the formulation (with changed parameters), touch **Start**.

Print example:

```
Component1
IdCm1
            39.2 g
Comp1 +
IdCm2
         Component2
Comp2 +
            23.3 g
  :
  :
  :
IdCm5
         Component5
            32.4 g
Comp5 +
               5
nDef
          117.6 g
Sum
```

Weighing in Percent


Purpose: This application is used to determine the percentage share or the percentage


difference of the sample related to a reference weight.

Options: You can enter a fixed reference weight or a reference percentage.

Combination options: Checkweighing, timer controlled functions, totalizing, statistics, formulation, second tare, individual identifiers, (SQmin function, DKD measurement uncertainty)

Configure Task: Weighing in Percent

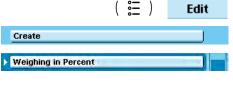


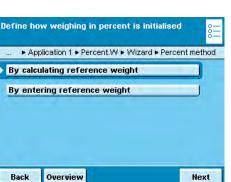

▶ Define how the weighing in percent will be initiated.

By calculating reference weight: The reference weight is determined by weighing a reference object.

By entering reference weight: The reference weight is entered as a numerical value.

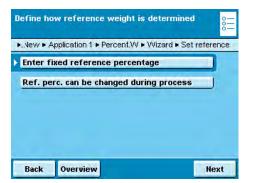


If the reference weight **should be** calculated.

▶ Define whether the reference percentage should be fixed or variable.

Enter fixed Reference percentage: You must now enter this.


*Reference percentage can be changed during process: You can enter this when carrying out the task.

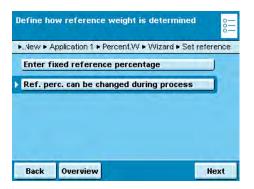

If the reference weight should be fixed:

Enter Fixed reference weight: You must enter this now.

*Reference weight can be changed during process: You can enter this when carrying out the task.

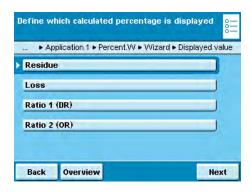


► Define which percent calculation should be used. **Example**: Weight value = 10 g, reference weight = 50 g

*Residue: The result is the percentage share of the weight value of the reference weight (Rest: Weight value/reference weight * 100%;in the example: 20%)


Loss: The result is the percentage loss between the weight value and the reference weight (loss: weight value - reference weight / reference weight * 100%; in the example: -80%)

Ratio 1(DR): The result is the percentage ratio of the difference to the weight value (Ratio 1: reference weight - weight value / weight value * 100%; in the example: 400%)

Ratio 2 (OR): The result is the percentage ratio of the reference weight to the weight value (Ratio 2: reference weight / weight value * 100%; in the example: 500%)

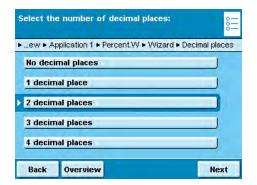

▶ Define the number of decimal places to be used for the display of the percentage. **none** ... **1**, ***2**, **3**, **4**

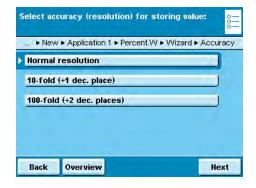

▶ Define the accuracy to be used to apply the weight value for the reference weight.

*Normal (display acc.) 10-fold (+1 dec. place) 100-fold (+2 dec. places)



- > An overview of all settings is displayed.
- Check all settings and change as required.
- ➤ A message asks whether you would like to add another application.
- ▶ If required, select additional applications to be combined.
- > Prompts for configuring weighing and printing appear.
- Check all settings and change as required.
- Enter a short name and a description for the new task.
- To save the new task, touch **Save**.

Execute Task: Weighing in Percent



- ▶ If you haven't already done so, go to Application Management.
- ➤ The task selection is displayed.
- ► Touch the desired task.

or

Start

If the desired task is already selected, touch **Start**.




- ▶ If required, enter the reference percentage (e. g. 100%) or the reference weight.
- ▶ Place the sample on the weighing pan.

In the example it is the reference sample.



- ➤ The weight value of the reference is displayed.
- ► To begin the Weighing in percent, touch **Start**.

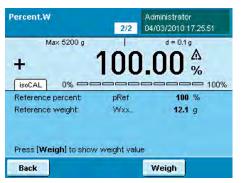

- ➤ The values for the reference percentage and the reference weight are displayed.
- ▶ Place the next sample on the weighing pan and touch **Weigh**.



- > The weight value is displayed.
- To toggle between the display of weight and percent, touch **Percent** or **Weigh**.

Timer-controlled Functions

Purpose: This application is used to trigger individual balance functions automatically at

a specific time or after specific time intervals. The time or interval must be within

24 hours (1 day).

Options: You can select the following functions:

Acoustic signalLock in readout

- Automatical printout of the display value

- Transfer of components for totalizing, formulation or statistics

Example: In order to determine the amount of evaporation of a sample over time, you can place the sample on the weighing pan and print out the weight value at fixed

intervals (e. g. every two minutes).

Combination options: Totalizing, statistics, formulation, mass unit conversion, second tare, memory,

individual identifiers (SQmin function, DKD measurement uncertainty)

(∷) Edit Create Timer-Controlled Funct. Define which function is triggered by the time New ▶ Application 1 ▶ Timer Ctr. ▶ Wizard ▶ Timer function Acoustic signal (beep tone) Lock in readout (hold display) Automatic printout of values Back Overview Next Please select the timer mode: New ▶ Application 1 ▶ Timer Ctr. ▶ Wizard ▶ Timer mode Timer set to intervals Timer set to a specific time Overview Next Back Please select the time input: . ▶ New ▶ Application 1 ▶ Timer Ctr. ▶ Wizard ▶ Time setting Enter fixed time Time can be changed during process

Overview

Configuring Task: Timer-controlled Functions

Define which function should be timer controlled.

*Acoustic signal (beep tone) The balance makes a beeping sound.

Lock in readout: The measured weight value

remains on the display for the defined time period.

Automatic printout of values: The displayed value is printed at the defined time.

Select the mode.

*Timer set to intervals: The function is triggered at regular intervals (e.g. every 30 sec.).

Timer set to a specific time: The function is triggered at the fixed time (after the start of the application) (e.g. at 08:00:00).

► Specify when the time and/or interval should be entered.

Define fixed time: You must enter this timepoint and/or interval now.


*Time setting can be changed: The time and/or interval is only entered when the task is carried out.



► Specify whether or not there should be an automatic restart after the timer-controlled function has been triggered.

*Off (donot restart)

On, restart automatically: The timer is restarted as soon as the timer controlled function has been carried out.

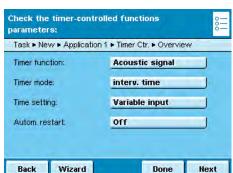

- ➤ An overview of all settings is displayed.
- Check all settings and change as required.
- ▶ A message asks whether you would like to add another application.
- ▶ If required, select additional applications to be combined.
- Now you will be prompted to configure the weighing and printing functions.
- Check all settings and change as required.
- Enter a short name and a description for the new task.
- To save the new task, touch Save.

Execute Task: Timer-controlled Functions

(등)

Start

- ▶ If you haven't already done so, go to Application Management.
- ➤ The task selection is displayed.
- ► Touch the desired task.

or

If the desired task is already selected, press Start.

If required, enter the time or interval.

(Enter hh:mm:ss, e. g. 00:00:30 for the 30-second interval)

► Follow the instructions on the display.



Back

Totalizing

Purpose: Adding up weight values and calculating values. The weight values of all individual

components are recorded and can be logged. The weight values and the component counter are saved in protected memory so that an interrupted totalizing process can be continued after the balance has been turned off or after a power failure.

Options: You can save a maximum of 65535 components.

User rights: All settings for the "Totalizing" application are saved under the active user profile.

Each user can make their own settings for this application. For this reason, make

sure that the desired user profile is selected before you start.

Results: When a user creates a local totalizing task, the results are only saved

locally for that user. The results of global tasks, however,


Combination options: Calculation, averaging, weighing in percent, parts counting, checkweighing,


are available to all users and the administrator.

mass unit conversion, second tare memory, individual identifiers, (SQmin function,

DKD measurement uncertainty)

(늘) Edit Create Totalizing

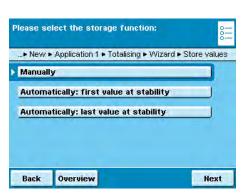

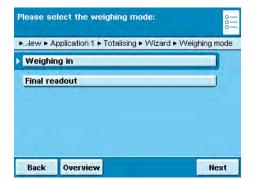

Configure Task: Totalizing

(* = factory setting)

▶ Define how the weight value will be applied.

*Manually: The weight value is applied as soon as the corresponding button is touched.

Automatically: first value at stability: The weight value is applied automatically as soon as the balance has stabilized after filling and the value exceeds the minimum load.

Automatically: last value at stability: The last weight value with stability is applied automatically, as soon as the balance is unloaded and the value falls short of the minimum load. If stability was reached repeatedly between measurements, the last stable weight value will be stored.

You must now select the minimum load to be applied:


(none, 10, 20, ...*100 .. 1000 digits).

Weighing Mode with Automatic Storage:

▶ Select the weighing mode to be used for storing the weight values.

Initial weighing: Initial weighing of samples in a container.

Finalreadout: Final weighing of samples in a container.

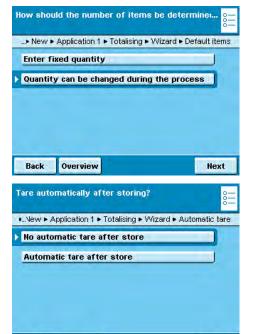


▶ Define when the number of measurements (items) will be defined.

Enter fixed quantity: The quantity of items is defined beforehand.

The quantity can be changed during the process

Enter the quantity of items when you execute the task.

▶ Define whether or not automatic taring should take place after storing the

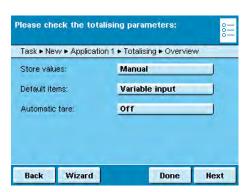

Back

Overview

- ➤ An overview of all settings is displayed.
 - Check all settings and change as required.
- > A message asks whether you would like to add another application.
- ▶ If required, select additional applications to be combined.
- Now you will be prompted to configure the weighing and printing functions.
- Check all settings and change as required.
- Enter a short name and a description for the new task.
- To save the new task, touch **Save**.

Execute Task: Totalizing

(📒)


Next

- ▶ If you haven't already done so, go to Task Management.
- ➤ The task selection is displayed.
- ► Touch the desired task

or

Start

- ▶ If the desired task is already selected, press Start.
- ▶ If required, enter the number of items that you want to add up.
- ▶ Place the first sample on the weighing pan.
- ► To start Totalizing, touch Start.
- During the task, you can show the results of previous measurements at any time by touching Calc.

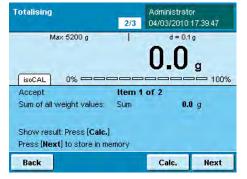

DKD Measurement Uncertainty

Purpose: This application is used to ensure that the measurement uncertainty is displayed

dynamically so that it conforms to the data documented on the DKD calibration

certificate.

Requirements: The function can only be used when the balance has been prepared for this before-


hand by a service technician. The service technician determines the measurement uncertainty of the balance via a DKD calibration onsite. The technician records on the DKD calibration certificate the measurements and the measurement uncertainty

for initial weighing. He also saves the calculated data on the balance.

Options: The measurement uncertainty can be displayed as an absolute value (U), a relative

value (U*) or as process accuracy (PA).

Combination options: Mass unit conversion, second tare memory, individual identifiers, (SQmin function)

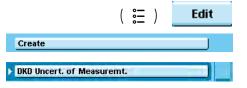

Configure Task: DKD Measurement Uncertainty

Select the measurement uncertainty display mode:


Select the measurement uncertainty display mode in relation to the maximum capacity.

Select the measurement uncertainty display mode in relation to the maximum capacity.

Absolute uncertainty: Absolute value (e. g. 12.00 g)

Relative uncertainty: Relative value (e. g. 0.000045%)

Process accuracy: e.g. 0.00013%

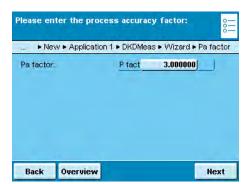


Next

Overview

Back

▶ If you have selected **process accuracy**, now enter the process accuracy factor.

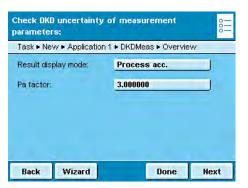

- An overview of all settings is displayed.
- ► Check all settings and change as required.
- ▶ A message asks whether you would like to add another application.
- ▶ If required, select additional applications to be combined.
- Now you will be prompted to configure the weighing and printing functions.
- Check all settings and change as required.
- Enter a short name and a description for the new task.
- To save the new task, touch Save.

Execute Task: DKD Measurement Uncertainty



- ▶ If you haven't already done so, go to Application Management.
- ➤ The task selection is displayed.

value set by a service technician.

Touch the desired task.

or

Start

- ▶ If the desired task is already selected, press **Start**.
- DKDMeas [S]: Administrator 09/03/2010 12:30:31 Max 5200 g d = 0.1g+ isoCAL U 0.08 g Cal./Adj. Menu DKDMeas [S]: Administrator 09/03/2010 12:31:14 Max 5200 g + isoCAL U* 1.4 % 0% ----Cal./Adj. Menu **DKDMeas** [S]: Administrator Max 5200 o
- ▶ If you have selected relative uncertainty the relative uncertainty of the

measurement is displayed in a tab as a percentage of the weight value.

If you have selected **Absolute uncertainty**, the uncertainty of the weight

value is displayed in a tab (e. g. U 12.01 g). This value is based on the DKD

▶ If you have selected **process accuracy**, the process accuracy is displayed in a tab in percent.

Cal./Adj.

isoCAL PA 4.3 %

Menu

0% ----- 100%

Second Tare Memory (Preset Tare)

Purpose: This application is used to define the second tare value. As soon as a second tare

value is used, **Net1** appears on the display for the respective net value.

Options: You can either use a weight value or enter a numerical value as a second tare value.

Combination options: Mass unit conversion, individual identifiers, (SQmin function, DKD measurement

uncertainty)

Configure Task: Second Tare Memory



- ▶ No pre-settings are required for this application.
- Now you will be to configure the weighing and printing functions.
- Check all settings and change as required.
- ▶ Enter a short name and a description for the new task.
- To save the new task, touch Save.

Execute Task: Second Tare Memory (Preset Tare)

(등)

- ▶ If you haven't already done so, go to Application Management.
- > The task selection is displayed.
- Touch the desired task.

or


Start

- If the desired task is already selected, touch Start.
- ▶ You have two options for defining the second tare.

Setting a Weight Value as Second Tare

▶ Place the object (container) on the balance whose weight value you want to use as the second tare.

Touch T1.

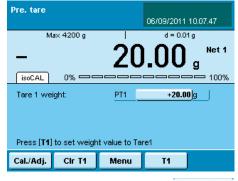

➤ The stored weight value is displayed as the second tare value. **Net 1** is also displayed next to the current weight value.

CIr T1


T1

Clr T1


To delete the tare memory, touch Delete T1

Entering a Numerical Value as a Second Tare

- ► Enter the tare value for **T1** (e.g. 20 g)
- ▶ Touch **T1** 0.00 g
 - ► Enter the desired value PT1 (e.g. 20)

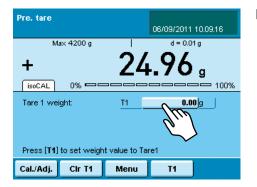

➤ The tare value is displayed.

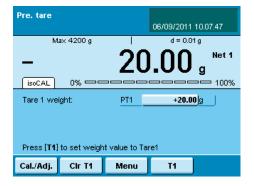

► To delete the tare value, touch **Delete T1**

Parts Counting

Purpose: This application is used to determine the number of objects which each have

approximately equal weight. This saves you from having to count

individual parts.


Options: You can calculate or enter the piece weight of individual objects.

Combination options: Checkweighing, timer-controlled functions, totalizing, statistics, formulation, mass

unit conversion, second tare memory, individual identifiers, (SQmin function, DKD

measurement uncertainty)

Configure Task: Parts Counting

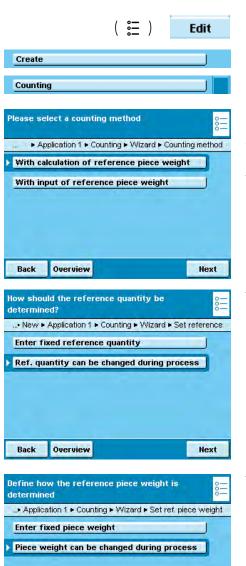

Overview

Next

► Select the method that you want to use.

*With calculation of piece weight: The piece weight is determined by weighing a known reference piece quantity.

Specification of piece weight: The piece weight is entered as a numerical value.

With calculation of the piece weight:

▶ Define how the reference piece count should be entered.

Enter fixed reference piece count: You must enter this reference piece count **nRef** now.


*Reference piece count can be changed during the process:

The reference piece count is entered when the task is executed.

When **entering** the piece count:

▶ Define how the piece weight should be entered.

Enter fixed piece weight: You must enter this piece weight **wRef** now. *Piece weight can be changed during process: The piece weight is entered when the task is carried out.

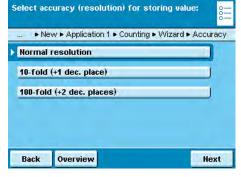

Define the accuracy to be used for the weight value.

*At normal resolution: As displayed

10-fold (+1dec. place): With one decimal place more than in the display **100-fold (+2 dec. places):** With two decimal places more than in the display




Only with calculation of piece weight:

▶ Define whether or not the piece weight should be updated when counting.

Optimisation of f

By pressing key, after prompt

*Automatic updating

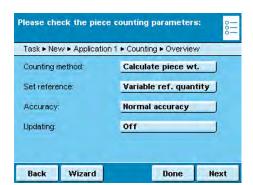


- ➤ An overview of all settings is displayed.
- Check all settings and change as required.
- ▶ A message asks whether you would like to add another application.
- ▶ If required, select additional applications to be combined.
- Now you will be prompted to configure the weighing and printing functions.
- Check all settings and change as required.
- ▶ Enter a short name and a description for the new task.
- To save the new task, touch **Save**.

Execute Task:

Parts Counting with Piece Weight Calculation



▶ If you haven't already done so, go to Application Management.

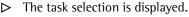

► Touch the desired task.

or

▶ If the desired task is already selected, touch **Start**.



- ➤ The last used reference piece quantity nRef is displayed.
- If you are working with a variable reference piece quantity, you must enter or change this now.
- ▶ Place this exact amount (nRef) of the objects to be counted on to the weighing pan.





- > The weight value for the reference objects is displayed.
- ► To determine the average piece weight, touch **Start**.



➤ The reference piece count and the average piece weight are displayed. You can now begin with the parts counting of unknown amounts.



- ▶ Place any amount of the objects on to the weighing pan.
- ➤ The determined piece count is displayed.
- To display the weight value of this amount, touch **Weigh**.



- ➤ The weighing results and all parameters are displayed.
- To count the next unknown amount, empty the weighing pan and place the next amount of objects on to it.
- > The calculated piece count and the weighing results are immediately updated.

Execute Task: Parts Counting by Entering the Piece Weight



- ▶ If you haven't already done so, go to Application Management.
- ➤ The task selection is displayed.
- ► Touch the desired task.


or

Start

If the desired task is already selected, touch Start.




- ➤ The last used piece weight wRef is displayed.
- ► Enter the piece weight of the reference object (e. g. 50 g).

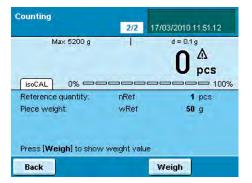

- ▶ The reference piece quantity and the entered piece weight are displayed.
- Place the sample with the objects to be counted on to the weighing pan.



- ➤ The determined piece count is displayed.
- To display the weight of the sample, touch **Weigh**.

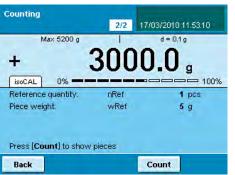

- ➤ The weight of the sample is displayed.
- To toggle between the display of weight value and piece count, touch **Count** or **Weigh**.
- To count the next quantity of objects, empty the weighing pan and place the next sample on to it.
- ➤ The weighing result is updated immediately.
- To display the piece count of this sample, touch **Count**.

Checkweighing

This application is used to check a weight value using preset control values. Purpose:


Options: Control values can be exact target values or tolerance range limit values within

which the check value must lie.

The check results are shown on the display. They can also be used for further

electronic editing by activating the control ports at the data output.

Totalizing, statistics, formulation, mass unit conversion, second tare memory, Combination options: individual identifiers, (SQmin function, DKD measurement uncertainty)

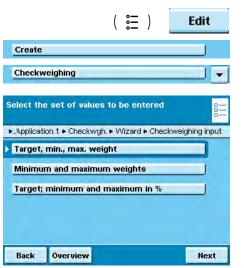


Configuring Task: Checkweighing

Define which values are to be used for checkweighing.

*Target, min., max. weight The target value and a tolerance range in absolute values

Minimum and maximum weights: Only a tolerance range in absolute values Target minimum and maximum in %: The target value and a tolerance range as a percentage

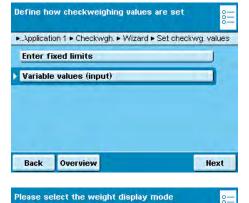

Absolute value

Back

Difference betw. target and actual

Overview

▶ Define how the control values will be entered.

Enter fixed limits: You must enter these values now.


*Variable values (input): The check values are entered when the task is carried

...New ▶ Application 1 ▶ Checkwgh. ▶ Wizard ▶ Display mode

Select the display mode for the results.

*Absolute value: Displays the absolute value.

Difference betw. target and actual: The deviation from the target value is then displayed.

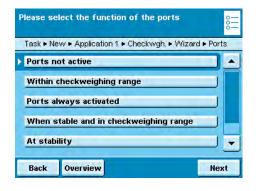

► Define how the ports should be activated depending on the check results *Ports not active


Within checkweighing range: Ports are only activated when the results are within the checkweighing range.

Ports always activated: The ports are always activated with each weight value. **When stable and in checkweighing range:** The ports are only activated as soon as the balance becomes stable and the results are in the checkweighing range. **At stability:** The ports are always activated as soon as the balance has stabilized. **Once at stability and in the checkweighing range:** The ports are activated on a one-time basis (not dynamically) as soon as the balance has stabilized and the results are in the checkweighing range; effective until the balance has been unloaded.


Define whether or not the automatic counter for positive results should be activated.

*No automatic counter

Counter for OK values: All positive checked results are counted.

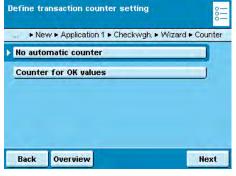

- An overview of all settings is displayed.Check all settings and change as required.
- ➤ A message asks whether you would like to add another application.
- ▶ If required, select additional applications to be combined.
- Now you will be prompted to configure the weighing and printing functions.
- Check all settings and change as required.
- ▶ Enter a short name and a description for the new task.
- To save the new task, touch **Save**.

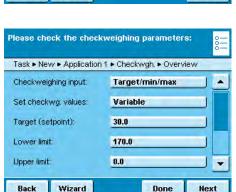

Execute Task: Checkweighing



- ▶ If you haven't already done so, go to Application Management.
- ➤ The task selection is displayed.
- Touch the desired task.

or

Start

If the desired task is already selected, touch **Start**.



Define check values by entry:

- ► Touch the input field Setp and enter the target.
- ▶ Touch the input field Min or Max and enter the minimum or maximum check values.

or

Моге

Define check values with current load:

- ▶ Place the weight on the weighing pan.
- Touch More.
- ► Touch the field of the check value whose current weight value you want to use.
- If required, repeat the process with additional weights for the other check values.
 To exit the input mode without saving the value, touch Back.

- To begin checkweighing, touch **Start**.
- > The entered check values are displayed graphically.
- Place the object to be checked on the weighing pan.



➤ The test results are displayed.

If the results are within tolerance, the weight results will always be displayed. In addition, they appear in green on in the graphic scale.



If the results are outside of the tolerance, the weight results are displayed in red (above maximum) or yellow (below minimum) on the graphic scale.



- ► To toggle the display to the check mode, touch **Check**
- Now the display will show whether the upper limit (HH) or the lower limit (LL) has been exceeded.
- ➤ To check additional objects, empty the weighing pan and place the next object on it.
- ▶ The display is updated dynamically.
- ► You can toggle between weighing and checkweighing at any time by touching **Weigh** or **Check**.





Importing/Exporting Data

You can use an SD memory card to transport and exchange data (import/export). This card is required when you want to save data externally or when you want to exchange data with other balances. For example, you can easily copy user profiles to several balances.

Inserting Memory Cards into the Display and Control Unit

The slot for the SD card is located at the back of the display and control unit.

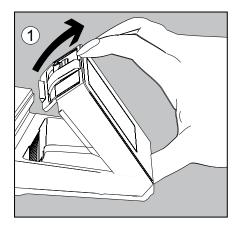

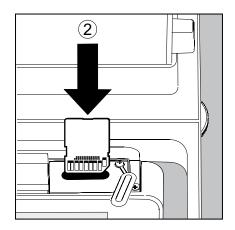

- 1. Tilt the control unit until is it nearly vertical.
- 2. Swivel the cover of the card slot to the front. Align the SD card so that the contacts are facing down.
- 3. Insert the card into the slot as far as it will go.
- ► To remove the card, press it against the resistance in the direction of the slot so that the card springs out.
- 4. Press the position retainer and swivel the display and operating unit back into the desired position.


Import/Export Data

The following data can be imported or exported via the SD card:

- Global tasks and separate tasks
- User profiles with their separate tasks
- Menu parameters
- Alibi memory
- Audit trail
- Calibration and adjustment data




You can open the data from the Alibi memory (xml Files) with any web browser and import them to Microsoft Excel as well.

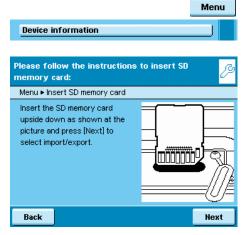


► Select the menu item Import/Export Data.

▶ Insert the SD card into the slot as shown in the picture.

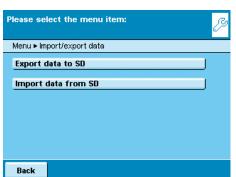

Select whether you wish to export or import data.

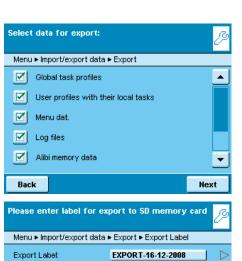

Export

► Select the data you wish to export and press **Next**.

- ▶ If required, change the name of the export folder and touch **Next**.
- ➤ The data is copied to the SD card.

An icon is shown on the display during the data transfer. The device cannot be operated during this time.


Back

Next

Import

In order for the data to be detected by the balance when importing them, you should not copy individual files, but always the entire subdirectory.

- data
- all
- EXPORT-dd-mm-yyyy
- task

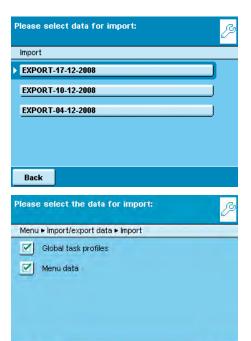

Back

► Select the folder on the SD card from which you want to import data.

- ▶ Select the data that you want to import and touch **Next**.
- ➤ The data will be imported from the SD card.
- ➤ An icon is shown on the display during the data transfer. The device cannot be operated during this time.

If the following error message appears:

"Wrong XML version. Please check the data."

Touch Next.

Next

- ➤ All previous data will be saved in the new format.
- ▶ Check your settings after importing the data (Menu, TASK and USER).
- ▶ If the XML files continue to cause problems, contact your local Sartorius service center.

Calibration and Adjustment

Background

During calibration, a check weight is used to determine how much the displayed value deviates from the actual measurement value. This deviation is compared with an entered target value and this deviation is then eliminated by a subsequent adjustment of the balance. During the Linearization, the deviation of the values from the ideal characteristic curve is cleared.



Not all functions/settings can be selected on balances used as legal measuring instruments!

When and how often

Calibration/Adjustment should be carried out on a regular basis, e. q. daily after the balance is turned on. In addition, it should be done after each leveling and always when the ambient conditions have changed (temperature, humidity or air pressure), and when the balance has been set up in another location.

isoCal

All Cubis balances are equipped with the automatic calibration function isoCal (time and temperature-controlled). You can adapt this function according to your requirements. You can set up a memory function that triggers a prompt for calibration/adjustment at specified intervals.




The process for the calibration/adjustment function may differ depending on the presets (see System Settings/Configuring Calibration/Adjustment). The factory settings have not been changed for the procedure described here.

Calibration/Adjustment Using Internal Check Weight

- If required, switch to operating mode.
- Make sure that the weighing pan is empty.
- Press TARE, to tare the balance.


(TARE)

Touch CAL.

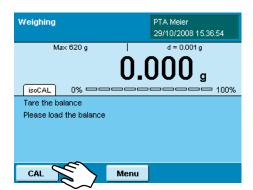



Touch Internal cal./adjustment and then Start.

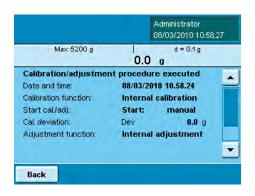

- The procedure will be executed.
- Wait until calibration/adjustment has been completed.
- > The results are displayed again after the procedure is completed.
- To return to operating mode, touch **Back**.

Calibration/Adjustment Using External Check Weight




(TARE)

An external check weight is required for this function. Please note the tolerance of the check weight being used.

- If required, switch to operating mode.
- ▶ Make sure that the weighing pan is empty.
- ▶ Press **TARE**, to tare the balance.
- ► Touch CAL



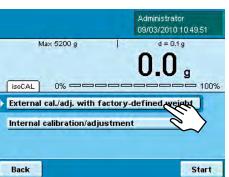

Calibration: Deviation

+ O.4 g


Calibration complete

Press [Exit] to exit without adjustment Press [Adjust] to activate adjustment

Exit Adjust

- Place the check weight on the balance.
- ► Touch Extternal cal./adj. with factory-defined weight.

- > Calibration is carried out (approx. 15 to 20 sec.).
- > After calibration, the determined deviation is displayed.
- ► To adjust the balance, touch **Adjust**
- To complete calibration without adjustment, touch Exit.

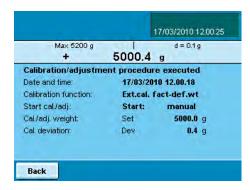

- ▶ Lastly, the results are shown on the display.
- ► To return to operating mode, touch **Back**.

User Management



You can configure the following settings in this menu:

- Creating new user profiles (only possible as administrator)
- Editing user profiles (modify, copy, delete, change and delete passwords, depending on user rights)
- Activate users

User Management

The user administration allows for defining of users with distinct sets of rights: the Administrator and multiple users.

The **Administrator** can use all functions (except those in Service) and he has all user rights. Only he can create new user profiles and assign individual rights. There is only one administrator.

A **user**, on the other hand, cannot use all functions. He has limited rights. Rights that are defined in the user profile. A max. of 15 user profiles can be created.

When the balance is first delivered, there are no user profiles, all settings can be carried out. The first user profile that is created is automatically an administrator. He can then create additional user profiles.



► To open the User Management menu, touch the USER (🖺)key.



> The menu opens and the selection of user profiles is displayed.


Creating User Profiles



This function is only available to an Administrator.

The following settings can be made in the user profile, they are only valid for that user:

- Language
- Display settings (color, brightness)
- Volume and acoustical signal on/off
- User rights
- Password protection
- User name and description.

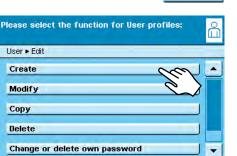

(음)

Edit

To create a new user profile, touch Edit.


► Touch Create.

The program now automatically guides you through the menu. Make your selections by touching the desired option, the program then takes you to the next prompt.

or

If the desired option is already selected (dark background), touch **Next**, to go to the next prompt.


If more than one option can be selected for a single setting, press **Next** after making your selection(s) to proceed to the next prompt.

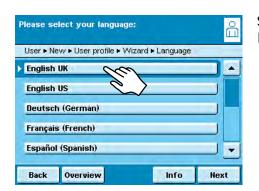

Back

Step 1: Select language

Touch the desired language.

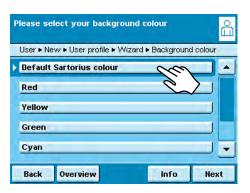

Steps 2 to 4: Setting the color and brightness of the display

- ► To select the background color, touch the desired color.

 The standard Sartorius factory setting for the background color is blue.
- ➤ The display changes to the color selection for the user field (see "Display in Operating Mode").
- ► Touch the desired color for the user field.

 The standard Sartorius factory setting for user fields is dark blue.
- ➤ The display changes to the brightness setting.
- ► To change the brightness of the display, touch the desired setting.

Step 5: Assign user rights

- ► Select the corresponding options to define which functions the user can access. The following rights can be assigned individually:
 - Creating, copying, modifying and deleting tasks
 - Importing data
 - Modifying the calibration and adjustment settings
 - Modifying the balance settings
 - Creating, modiying and deleting timer-controlled actions
 Administrators can access all functions.
- ► To move to the next menu item, press **Next**.



Step 6: Define password protection


If you assign a password, this profile cannot be used until the password has been entered.

Select the corresponding option to define whether the user profile is password-protected or not.

The password will be defined in the step after next.


Only the Administrator can open the user profile without the password (for example, if the user forgets the password) and assign a different password. The Administrator can remove the password protection so that a new password can be set, if desired.

If the administrator forgets the password: Please contact your Sartorius Service Center

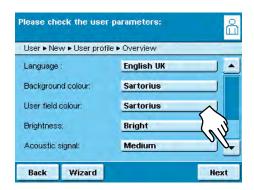


Step 7: Check all settings

- ▶ Use the scroll bar or press the down arrow to view all settings.
- ► To modify a setting, press the corresponding button. This opens a screen showing the options for that setting.
- ▶ If the settings are correct, press Next.

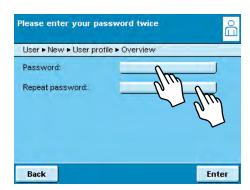


Step 8: Define the password (If password protection enabled)

- ▶ If you selected the option for password protection in Step 7 (see above), you are now prompted to enter the desired password twice.
- Press the upper button.
- ► Enter the desired password (minimum 4 characters, maximum 8; only numbers and uppercase letters are permitted).
- Press the lower button, and enter the password again.
- Touch Next.

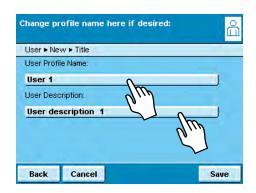

Step 9: Enter user names and descriptions

- To enter a name for the user profile, touch the **User name** button. The user name is shown on the top right of the display during operation as long as the respective user profile is active.
- ► Enter the name (max. 40 characters) and touch **OK**.
- ➤ To enter a description of the user profile, touch the **User Description** button. The user description can be viewed at any time during operation of the balance by pressing the user info area in the display (see "User Interface (Touch Screen)").
- ► Enter the description and touch **OK**.
- ► To complete the process, touch **Save**.

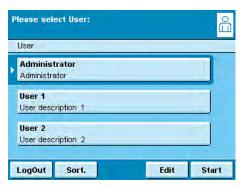

- ➤ The program switches to user selection. The newly created user profile is placed at the end of the list.
- ▶ When you want to activate a new user, just touch the button with his name.
- ➤ The program switches to the operating mode and the selected user is activated.

or

If you want to set up additional user profiles, touch Edit. and repeat the steps described.

or

▶ When you want to switch to operating mode without switching the user, touch **Start**.

Activating Users




With the balance in the operating mode, the active user is displayed in the user field at the top right of the screen.

To activate another user, you must switch to the **User Management** menu.



ightharpoonup Press the USER key ($\stackrel{\circ}{\square}$).

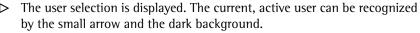

- Touch the user you want to activate.
- ▶ If the user is password-protected, enter the password and confirm the entry.
- > The program switches back to operating mode with the selected user activated and displayed at the top right of the user field.




Editing User Profiles




6

- ► If you have not already done so, open the User Management menu by touching the USER () key.
- To go to the Edit menu, touch **Edit** in the user selection.
- > The selection of edit functions is displayed.

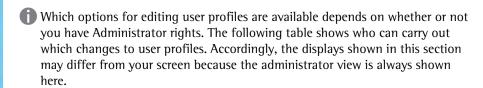

Option	Administrator	User
Change own profile	\sqsubseteq	\square
Change other profile	\sqsubseteq	
Copy profile	\sqsubseteq	
Delete profile	\sqsubseteq	
Change or delete own password	\sqsubseteq	\square
Change or delete other user passwords	\sqsubseteq	

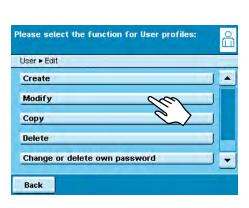

Please select User:

Administrator Administrator

User description 1


User

User 1

LogOut


Changing User Profiles

- Select Modify.
- If you are an administrator, you must now select the profile that you want to change and touch **Next**.
- > The overview of all settings for this profile is displayed. You can now select and change the parameters one by one.
- If you want to use the automatic user guide (step-by-step), touch **Wizard**. Now all setting options will be displayed one by one to be changed.
- Touch Next.
- ▶ If required, change the user name and description.
- Touch Save.
- ➤ The user profile has been changed. The display changes to the user selection.

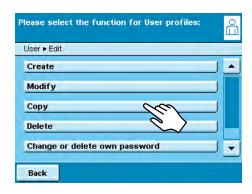

Copying User Profiles

- This function can only be carried out by an administrator.
- When you copy a user profile with password protection, the password is not transferred to the copy, rather it is deleted.
- Select Copy.
- ➤ All available user profiles are displayed.
- ► Select the profile that you want to copy.
- Touch Next.
- ▶ Enter a new name and description for the copied user profile.
- Touch Save.
- ▶ The display switches to the user selection in which the new profile is displayed.

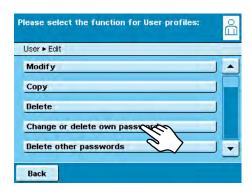

Deleting User Profiles


- This function can only be carried out by an administrator.
- ► Select **Delete**.
- ➤ All available user profiles are displayed.
- Select the profiles that you want to delete and touch Next.

or

- If you want to delete all profiles, touch AII.
- ➤ All profiles selected for deletion are displayed in a security prompt.
- ▶ If you want to change your selection, touch **No.**

or

Confirm the deletion with Yes.

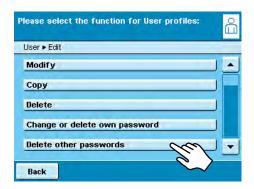


Changing or Deleting Own Password

- ▶ If you haven't already done so, activate your own user profile.
- ► Touch **Edit.** in the user selection.
- Select Change or delete own password.
- ► To change the password, touch Yes.
- ▶ Touch both buttons one after the other and enter the new password twice.

or

- ► To **delete** the password, touch **No**.
- ► Touch Next.
- ➤ The display switches to the selection of change functions.

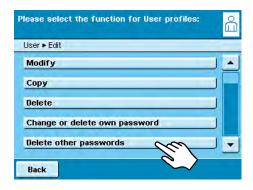

Deleting Other User Passwords

This function is only available to an Administrator.


- Activate your own profile.
- ► Touch **Edit.** in the user selection.
- Select Change or delete other password.
- ➤ An overview of all users with passwords is displayed.
- Select all users whose passwords you want to delete and touch Next.

or

- If you want to delete all passwords, touch AII.
- ▶ All users whose passwords you selected are displayed in a security prompt.
- If you want to change your selection, touch No.
- Confirm the deletion with Yes.
- ➤ The display switches to the overview of user settings.

Interfaces

Purpose

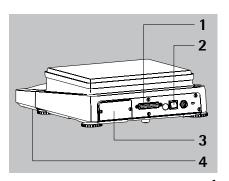

Interfaces are used to exchange data with ported peripheral devices:

Measured values and calculated values can be output to a printer, PC or control display; conversely, control commands and data inputs can be sent to ported devices (PC, keyboard, foot switch, bar code scanner).

Except for PS2, each interface has to be configured according to the peripheral device and desired function. No error messages are generated just because no device is connected to an interface port (open data port).

Cubis series balances have at least three interfaces:

Peripheral port (25-pin interface), Com A

- 2 USB interface for a PC connection, Com B
- 3 The slot may contain other ports (Com C):
- 25-pin interface (YD001MS-R)
- 9-pin interface with 6-pin PS2 port (YD001MS-P)
- Bluetooth® Module (YD001MS-B)
- 4 Ethernet interface (on the bottom of the display unit), Com D

Data Exchange Protocols

For data exchange, the interfaces are configured with the following protocols:

- Printer output
- SBI (Sartorius Balance Interface): Sartorius standard protocol for output to a PC or control unit. This simple ASCII-based protocol allows you to use ESC commands from your PC to control the basic weighing functions.
- xBPI (eXtended Balance Processor Interface, also called X-Bus): binary protocol
 with extended command volume. This protocol lets you control numerous
 weighing functions. For further information on this, please contact Sartorius.
- SICS (Standard Interface Common Set): This interface protocol enables operation and control of the balance via a connected PC. You can read out measurement data, enter weighing commands and activate all operating functions.

To use the protocols, application software must be installed on the PC, e.g. SartoTerminal.

Synchronization

During data communication between balance and PC, messages consisting of ASCII or binary characters are transmitted via the interface. For error-free data exchange, parameters for baud rate, parity, handshake mode and character format must be identical for both units.

You can configure the respective settings in the System Settings menu. In addition to these settings, data output for the balance can also be made dependent on several conditions that are defined in the individual tasks. These conditions are described under each of the tasks.

USB Port (PC)

Purpose

Any Cubis balance can be connected to a PC equipped with a USB port. A virtual serial interface (virtual COM port) is set up as a device type at the USB port. This virtual serial interface is identified und operated by the application program. The protocols xBPI, SBI and SICS can be transferred via the USB port.



The USB port is designed for the laboratory environment and is not suitable for use in purely industrial environments. Full IP protection is only guaranteed when the USB cover is closed.

System Requirements

- Computer (PC) with Windows 98SE®, Windows ME®, Windows 2000®,
 Windows XP®, Windows Vista® or Windows 7®
- Available USB port on the PC
- USB cable

Software Driver and Installation Guides

The VCP driver, used to set up the virtual interface on the computer, can be downloaded from the internet: http://www.ftdichip.com/FTDrivers.htm

The installation guides for the drivers can be found here: http://www.ftdichip.com/Documents/InstallGuides.htm

Connecting the Balance via USB



The current USB port for the computer is established when the software driver is being installed. The driver must be re-installed every time you wish to change the port. Therefore, choose one USB port that can permanently or regularly be used to connect the balance.

- Switch off the balance.
- Unplug the balance from the mains.
- Connect the USB cable to the balance and to the USB port on the computer.
- Plug the balance into the mains again and switch it on.
- ▶ Windows detects the device connected to the USB port. If the device is being connected for the first time, the Windows Installation Wizard will run.

Installing Software Drivers

- Run the Installation Wizard for the driver.
- Follow the instructions that appear.
- To complete the installation, click on Finish.
- ➤ The virtual interface is now ready for operation.

Windows® usually adds the virtual port in the position following your highest-numbered COM port.

Example:

For a PC with up to 4 COM ports, the new virtual port would then be COM5 (see Device Manager).

Installation guides for Windows XP®, Windows Vista® and Windows 7®

Changing the Port No.

If you use the USB interface with a program that limits the number of COM port designations (e. g. only COM1, 2, 3, 4), you may have to assign one of these port numbers to the new virtual port.

- Open the setting for the **USB serial port** in the Windows® system control panel:
 - START > My Computer > Control Panel
 - System > Hardware > Device Manager
- Open the **Connections** submenu.
- Double-click on USB Serial Port.
- Select Port Settings / Advanced.

Change Latency Time

- Open the settings for the USB serial port, following the above instructions.
- For a faster rate of communication, change the setting for the **Latency Timer** to 1 msec.

Plug & Play Mode in Autoprint (SBI)

- Open the settings for the USB serial port, following the above instructions.
- Stop the **Plug & Play mode** from running.

Uninstalling the Driver

The software driver for the USB connection can be uninstalled with the Windows® Uninstaller.

PS2 Interfaces for Bar Code Scanner or Keyboard

Via the PS2 port you can connect a bar code scanner or a PC keyboard to enter data and operate the balance.

Installation Connecting the Device The interface installation is described in the installation instructions supplied.

- Plug the connector plug into the PS2 port on the balance.
- ➤ The input device is operational, no further settings are required.

Using the Input Device

You can input entries directly via the keyboard or bar code scanner whenever an input field is open in the display of the balance.

Menu Calc. Start F3 F4 F5

Param.

F2

CAL

F1

Key Assignment of the PC Keyboard

The five context-dependent buttons located at the bottom edge of the display are operated via the Function keys F1 to F5 on the PC keyboard.

Back Esc

Whenever the last button on the right is assigned with the function **Back**, this function can also be activated using the Esckey on the keyboard.

F9

The **TASK** key is operated via **F9** (Open Task menu).

F10

The **USER** key is operated via **F10** (Open User menu).

Print


The **PRINT** key is operated via the Print key of the keyboard.

Shift F1

Use the shift F1 to select the **first entry** of a menu.

Shift F2

Use shift F2 to select the **second entry** of a menu. etc.

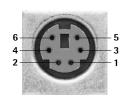

Pin Assignment PS2

Pin 1: Keyboard_DATA (universal)


Pin 2: Not assigned (larger)

Pin 3: Internal ground (GND)

Pin 4: +5 V


Pin 5: Keyboard_CLK (Set)

Pin 6: Not assigned (smaller)

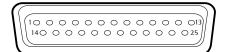

Interfaces (RS-232) 25-pin and 9-pin

The balance is standard-equipped with a 25-pin interface (**peripheral port**) where different Sartorius peripherial devices can be connected, e. g. hand or foot switch. This interface is also used for the control lines for the **Checkweighing** application.

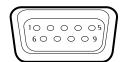

A 9-pin interface can be optionally installed. This port is intended for connecting a PC.



Warning When Using Pre-wired RS-232 Connecting Cables for the 25-pin Interface:

RS-232 cables purchased from other manufacturers often have incorrect pin assignments for use with Sartorius devices. Failure to do so may damage or destroy your weighing system and/or peripheral devices!

- ▶ Be sure to check the pin assignments before you connect cables purchased from other manufacturers.
- Disconnect any lines assigned differently (e. g. Pin 6).

Features (*= Factory settings)

Type of interface:	Serial Interface		
Interface operating mode:	Full duplex		
Level:	RS-232		
Interface connector:	D-SUB female connector, 25- or 9-contact		
Transmission rate:	600, 1200, 2400, 4800, *9600 and		
	19200 baud (selectable)		
Parity:	*Even, odd, none; blank spaces (selectable)		
Character transmission:	Start bit, 7/8* bit ASCII, parity, *1 or 2 stop bits		
	(selectable)		
Handshake (selectable): 25-pin	For 2-wire interface: Software (XON/XOFF)		
•	for 4-wire interface: *Hardware (CTS/DTR)		
9-pin	Hardware (CTS/RTS)		
Operating mode:	*SBI, xBPI, printer, SICS, 2nd display		
Manual print mode	Without stability, *after stability		
Auto print mode	*Without stability, at stability, after load change		
Cancel automatic printing:	Not possible		
Time-dependent automatic			
printout:	After 1 display update		
Data output of balance:	16 or *22 characters		
Tare after individual printout:	Off		
Basic values, application:	Off		

Connectable Peripherials

The following devices can be ported to the interface:

- Verifiable data printer YDP03-0CE, YDP10-0CE, YDP20-0CE
- Verifiable data printer with Bluetooth data transmission YDP10BT-0CE (module required)
- Universal (serial printer)
- Strip/label printer YDP04IS-OCEUV
- Hand switch YHS01
- Foot switch YFS01, YPE01RC
- External control display YRD11Z
- Remote display YRD03Z


Pin Assignment Chart for Connecting Peripherals

For Sartorius peripheral devices Purpose:

Female Interface Connector: 25-pin D-Submini (DB25S) with screwed fastening

Required Male Connector (Recommended): 25-pin D-Submini, DB25S, with integrated shielded cable clamp and shield plate

assembly (Amp type 826 985-1C) and fastening screws (Amp type 164 868-1)

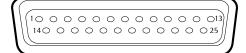

Pin assignment 25-pin socket, RS 232:

Pin 1: Signal ground Pin 14: Internal ground (GND) Pin 2: Data output (TxD) Pin 15: Control input/output 12) Pin 3: Data input (RxD) Pin 16: Control input/output 2²) Pin 4: Signal GND Pin 17: Control input/output 32) Pin 5: Clear to Send (CTS) Pin 18: Control input/output 42) Pin 6: Not used Pin 19: Control input/output 5²) Pin 7: Internal ground (GND) Pin 20: Data Terminal Ready (DTR)

Pin 8: Internal ground (GND) Pin 21: Not used Pin 9: Not used Pin 22: Not used Pin 10: Not used Pin 23: Not used Pin 11: + 12 V output Pin 24: Not used Pin 12: Reset _ Out1) Pin 25: +5 V output

Pin 13: + 5 V output

= Peripherals restart

²) = Allocation of control input/outputs menu-configurable (see below)

Control Inputs/Outputs

Data Output / Setting	Pin 15	Pin 16	Pin 17	Pin 18	Pin 19
Infrared sensor YHS01MS	Input 1: *"PRINT" key	Output 2: ** "Smaller"	Output 3: ** "Equal"	Output 4: ** "Larger"	Output 5: **"Setp"
Hand switch YHS02	Input 1: *"PRINT" key	Output 2: ** "Smaller"	Output 3: ** "Equal"	Output 4: ** "Larger"	Output 5: **"Setp"
Foot switch YFS01	Input 1: *"PRINT" key	Output 2: ** "Smaller"	Output 3: ** "Equal"	Output 4: ** "Larger"	Output 5: **"Setp"
Triple Foot switch YPE01RC	lnput 1: Left *"PRINT" key	Input 2: Right *Left Shieldpl.key	No function (disabled)	No function (disabled)	Input 5: Center *"TARE" key
Control display YRD11Z	Output 1: Currently no function	Output 2: ** "Smaller"	Output 3: ** "Equal"	Output 4: ** "Larger"	Output 5: **"Setp"
Control inputs	lnput 1: *"PRINT" key	Input 2: *Left Shieldpl.key	Input 3: *no function	Input 4: *no function	Input 5: *"TARE" key
Control outputs	Output 1: Currently no function	Output 2: ** "Smaller"	Output 3: ** "Equal"	Output 4: ** "Larger"	Output 5: **"Setp"

Default allocation of input, otherwise configurable

137


Allocation of outputs, on checkweigher

Cabling Diagram 25-pin Interface


Diagram for connecting a computer or other peripheral device to the balance using the RS232/V24 standard and cables up to 15 m (50 ft.) long



Do not connect any other pins to the cable connector of the balance!

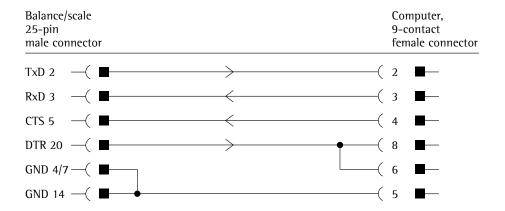

Cable type: AWG 2 specification

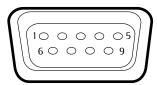

Pin assignment 9-pin socket, RS-232:

Pin 1: Not used

Pin 2: Data output (TxD)

Pin 3: Data input (RxD)

Pin 4: Not used

Pin 5: Internal ground (GND)

Pin 6: Not used

Pin 7: Clear to Send (CTS)


Pin 8: Request to Send (RTS)

Pin 9: Not used

Establish a connection via a conventional RS-232 cable.

Configuring Serial Ports



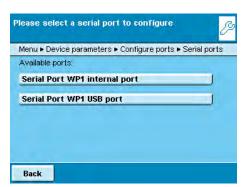

- The available interfaces are displayed.
- Select the interface you want to configure.
- Determine all settings for this interface.
- To save settings, touch **Save**.

Configuration Options for the Serial Interfaces

The following **operating modes** are possible for these interfaces:


- No function; the interfaces is blocked
- xBPl
- SB1
- Remote display (SBI)
- Printer output
- SICS



Factory settings are marked with an *.

Configuration Options for the xBPI Operating Mode


For this operating mode, you can only configure baud rate and number of stop bits.

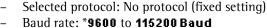

Data bits: 8 (fixed setting)

Parity: Odd (fixed setting)

Stop bits: *1 or 2 stop bits

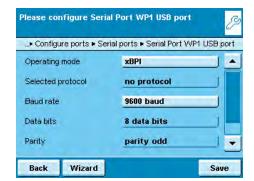

Configuration Options for the SBI Operating Mode and Remote Display

You can change all settings for this operating mode.


-▶ Configure ports ▶ Serial ports ▶ Serial Port WP1 USB port Selected protocol: No protocol/software handshake XON;XOFF/*Hardware handshake RTS;CTS
 - Baud rate: **300** to **19200 baud** (for remote display up to 115200 baud), *9600
 - Data bits: 7 or *8
 - Parity: None/*Odd (odd)/Euen (even)
 - Stop bits: *1 or 2 stop bits Log data: *Turn of f/Turn on




SBI

Please configure Serial Port WP1 USB port

Operating mode

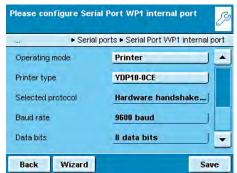

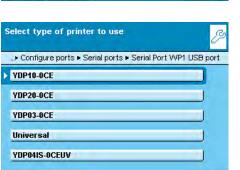

Back

OK


Configuration Options for the Printer Output Operating Mode

You can modify all settings for this operating mode depending on the configured printer type.

- Printer type: YDP10-0CE/YDP20-0CE/YDP03-0CE/Universal/ YDP04IS-0CEUV
- Selected protocol: Hardware handshakeBaud rate: 300 to 19200 baud (*9600)
- Data bits: 7 or *8
- Parity: None/*Odd (odd)/odd (even)
- Stop bits: *1 or 2 stop bits
- Log data: *Turn of f/Turn on

The following setting options are available for the **Printer types**:

- YDP10-0CE: No setting option, all parameters are permanently defined
- **YDP20-0CE**: All can be configured except stop bits
- YDP03-0CE: All can be configured except data bits and stop bits
- Universal: All can be configured
- YDP04IS-0CEUV: No setting option, all parameters are permanently defined
- YDP10BT-0CE (only for optional Bluetooth module, Com C): No setting option, all parameters are permanently defined

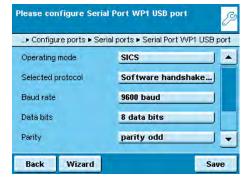

Configuration Options for the SICS Operating Mode

You can change all settings for this operating mode.

- Selected protocol: No protocol/*Software handshake XON;XOFF/*Hardware handshake RTS;CTS
- Baud rate: 300 to 115200 Baud, *9600
- Data bits: 7 or *8
- Parity: None/*Odd/Even
- Stop bits: *1 or 2 stop bits
- Log data: *Turn of f/Turn on

Explanations for the Setting Parameters

Selected Protocol/Handshake

The SBI (Sartorius Balance Interface) has transmit and receive buffers. Data transmission can be operated with a hardware or software handshake

Hardware Handshake (CTS/DTR):

With a hardware handshake (4-wire interface) protocol, the handshake is controlled via the CTS and DTR lines.

Software Handshake (XON, XOFF) The software

The software handshake is controlled via the XON and XOFF control commands. When the device is switched on, XON must be transmitted to enable any connected device to communicate.

Data transmission:


Balance --- byte ---> --- byte ---> (receiving device) (transmitting device) --- byte ---> --- byte ---> <--- XOFF ------ byte ---> --- byte ---> (Pause) <--- XON ------ byte ---> --- byte ---> --- byte ---> --- byte --->

Transmitting Device: Once XOFF has been received, it prevents further transmission of characters.

Once XON has been received it re-enables the transmitting device to send data.

Receiving Device: To prevent too many control commands from being received at one time, XON

is not transmitted until the buffer is almost empty.

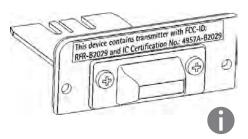

Bluetooth® Interface (COM C, Optional)


How to assemble the Bluetooth® module is described in the installation instructions supplied.

For transmitting data via the Bluetooth module, the data output has to be configured in the System Settings as follows.

Device information Configure data output Configure ports


Configuring the Bluetooth® Interface

- ▶ If appropriate, toggle to the System Settings menu.
- ▶ Select the **Device information** submenu and slide the right scroll bar down.
- Select Configure ports.
- Open the Serial ports submenu.
- Select Available ports.

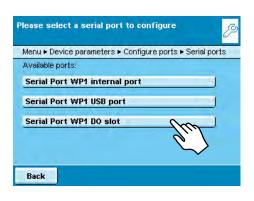


The following **operating modes** are available for this port:


- No function
- SBI (data exchange with PC or notebook)
- Printer output
- SBI (data exchange with PC or notebook)
- Select the desired operating mode

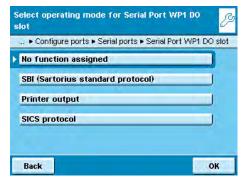

Bluetooth Printer Output Operating Mode

- ► If required, you can change the device name (only with Bluetooth modules from Version 04-10-01).
- To connect a PC/Notebook, select **PC or other**.

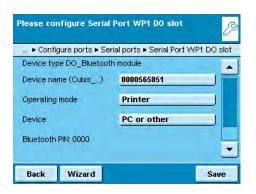

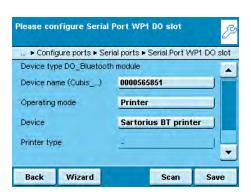

Connecting a Bluetooth Printer

- ▶ Make sure that the printer to which you wish to connect is ready to operate.
- ► Select the **Sartorius BT printer** as the device.
- To establish the connection, select **Scan**.
- ➤ The balance will now search for the Bluetooth device. This process may take time. Wait for the next message to appear on the display.



- ➤ A message will appear to confirm whether the printer has been found.
- ► To save settings, touch **Save**.
- Saving settings: Touch 2x on Back.
- ➤ The connection will initially be terminated after saving. The LED on the printer will continue to flash blue/red.
- SAs soon as the configuration menu is exited, the connection will be re-established. The LED on the printer will then illuminate permanently (blue).

Network Interface (Ethernet)

Purpose

This interface lets you integrate your balance into a TCP/IP-based network.

System Requirements

The person configuring the Ethernet interface should have basic knowledge of TCP/IP-base networks and network technologies in general.

The Ethernet interface is connected to the local network via the RJ45 slot.


Features

Transfer rate:	10 Mbs/sec (10BASE-T, Ethernet) and 100 Mbs/sec (100BASE-TX, Fast Ethernet), auto-recognition (10/100, HalfDX/FullDX)
Connection type:	Network protocol: ModBus/TCP
Transport:	TCP/IP or UDP/IP
HMS certificates:	Compatible with ModBus TCP standard of the ModBus Organisation, industry-compatible CE, UL, cUL
Cable:	Twisted pair, shielded, CAT-5 or higher, 1:1, UTP / STP, male connector RJ45; e. g. patch cable CAT5 depending on application (straight/crossover)
Cable impedance:	150 Ohm
Cable length to HUB:	Max. 30 m

Connecting Network Cables


The slot for the network port is located at the back of the control unit.


► Tilt the control unit (1) until it is nearly vertical.

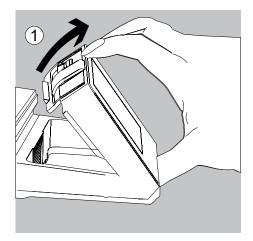


On the bottom panel of the display unit:

► Turn the locking lever (2) 90°.

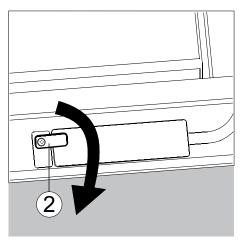


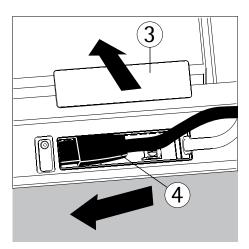

- Remove the cover plate (3).
- ▶ Insert the Ethernet cable plug (4) so that it clicks audibly into place.

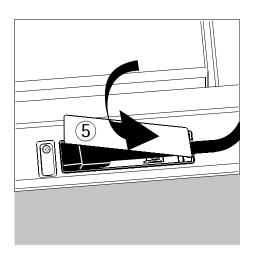

- Reattach the cover plate (5).
- ► Turn the locking lever back to its locked position.

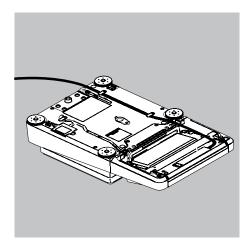



Exercise caution to avoid glass breakage for models with a draft shield.

- ► Turn over the balance and place it on a soft surface.
- ▶ Lay the Ethernet cable (6) through the cable channel of the weighing module.

Configuring Network Operation (Ethernet)

Preparation:

Before you can start with the configuration, you will require some information about your network. Please contact your network administrator. The important question is whether your network supports DHCP or not.

DHCP

With DHCP, you can automatically link a (new) balance to an existing network without having to configure it manually. Normally, the balance (client) only requires the automatic acquisition setting for IP address (factory setting). In this case, no additional information is required.


DHCP is supported DHCP is not supported

In this case, you need the following information from the network administrator:

- IP address
- Subnet mask
- Default gateway address
 Device name (host name): This name can be changed if required.
 This name is used to select the balance in the network.

wSelect Network Settings (Fixed Setting)


- ▶ If appropriate, toggle to the System Settings menu.
- ▶ Select the **Device information** submenu and slide the right scroll bar down.
- Select Configure ports.
- Open the Network parameters (Ethernet) submenu.

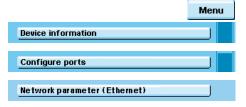

► To change the IP settings, touch the input field behind **Network configuration**input field.




Select Use fixed IP address and confirm with OK.

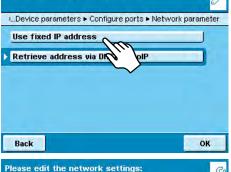

In the overview, check all settings and change as required.

Device name (host name): You can select this name freely. This name is used to select the balance in the network.

- ▶ If you then want to produce an Ethernet connection, write down the IP address now.
- To save the network configuration, touch **Save**.
- ➤ The Configure ports submenu is displayed.

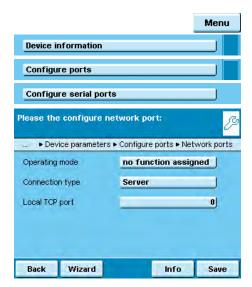

Configuring Network Interface (Com D)

- ▶ If appropriate, toggle to the System Settings menu.
- ▶ Select the **Device information** submenu and slide the right scroll bar down.
- Select Configure ports.
- ▶ Open the **Network ports**. submenu.

▶ Define the operating mode and, depending on the connection type, either Server port and address or the local TCP port.

Operating mode: No function/XBPI/SBI/Printer output/SICS protocol Connection type (of the balance): Server/Client

► To save settings, touch **Save**.

Examples:

For the connection type **Balance = Server** you only need to enter the local TCP port (e. g. for working with the Sartorius **SartoCollect** software for data communication)

For the connection type **Balance = Client**, you have to enter the server TCP port and the server IP address, (e. g. for working with "**Virtual serial port emulator"** as the server for all previous serial PC applications with the balance or for printing on a network printer server.)

Remote Control of the Balance from a PC via Ethernet

System Requirements

- PC with Internet browser
- Operating systems Windows, Mac OS, Linux, Unix and Solaris.

Installation

- Connect the balance to your network via the Ethernet interface.
- ► Enter the IP address of the balance in the address field of your Windows browser. The IP address of your balance is listed under Menu/Device Information/Basic data.

Port Release

- If you are having problems with the connection, you must configure your firewall to release the right TCP/IP UDP port.
- If you are using a router and want to access your computer via the Internet via your IP address, the port must also be set in the router's virtual server settings.
 This means that you must forward the port to the IP address of the server.
- If you still have problems, you may have to deactivate your firewall temporarily while using a browser connection.



Selecting Functions

- ► Click on the desired function in the browser.
- ► To exit a function, click on >> Back or on the Go back key of your browser.

Remote configuration (VNC): Use this function for remote control of the balance. The display screen on the balance is reproduced in the browser, you can operate the balance now by using the mouse and keyboard of your PC.

Software version: You can view the version of the balance, control unit and application software.

Screenshot: You can print the current screen or save it as GIF file.

Browse alibi memory: You can view the Alibi memory.

Show settings: You can view the following settings for log files:

- Task profiles
- User profiles
- System settings (menu)
- Service information (service)
- Interface information (ports)
- Timer controlled actions

Show log files: You can view respective log files.


Viewing a print file directly in an Internet Browser:

Prerequisites:

The configuration settings can be found in the chapter System Settings (Menu): "Print to File." Select the following file format for viewing directly in an Internet browser: Print to TXT file.

- In the browser, click on "Show Log Files" and then "Print to File."
- To view the printout directly in the Internet browser, click on "print.txt."
- ➤ The printout is displayed in the Internet browser.



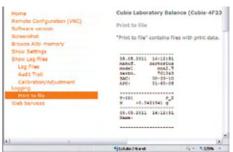

Data Output

You can define the data output parameter so that output is activated either when a print command is received or automatically synchronized with the display or at defined intervals (see application programs and autoprint settings).

Data Output by Print Command

The print command can be transmitted by pressing (\square) or by a software command (Esc P).

Automatic Data Output

In the **Autoprint** mode, data is output to the data interface port without an extra print command. You can have synchronized data output automatically at defined display update intervals, with or without the stability parameter. The interval time depends on the balance operating status and balance type.

If the automatic data output is activated in the Device Configuration, it starts immediately after the balance is turned on. You can also configure whether the automatic data output can be stopped and started by pressing the () key.

Data Output Formats

You can output the value displayed in the measured value line and the weight unit, with or without a data ID code. Configure this output parameter in the Device Configuration menu (Menu/Device Configuration/Configure data output/Line format).

Example: Output Without an ID Code

+ 253 pcs

16 characters are output

Example: Output Without an ID Code

Qnt + 253 pcs

22 characters are output

Data Output Format with 16 Characters

Display segments that are not activated are output as spaces. Display values without a decimal point are output without a decimal point. The type of character that can be output depends on the character's position.

Normal Operation

	- 1															
Position	1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16
	+	Α	Α	Α	Α	Α	Α	Α	A	Α	*	Е	Е	Е	CR	LF
or	-							•		•		*	*	*		
or		*	*	*	*	*	*	*	*	*	*					

*: Spaces

CR: Carriage return

A: Displayed characters

LF: Line feed

E: Characters for the unit

_				
۱n	ecia	1 ()	utn	uits

or*

Position	1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16
	*	*	*	*	*	*	_	_	*	*	*	*	*	*	CR	LF
or*		*	*	*	*	*	Α	*	*	*	*	*	*	*	CR	LF
or*		*	*	*	*	*	Α	В	*	*	*	*	*	*	CR	LF
and only	upc	n re	ques	t wit	h ES	C wo	(no	prin	t con	nmar	nd):					
		*	*	*	*	*	W	*	χ	Χ	Χ	Υ	Υ	Υ	CR	LF

*: Spaces A = C: Adjustment

AB = -: Final readout W: Draft shield status (optional):

X X

Χ

Manual draft shield operation

CR LF

A = H: Overload 1: lonizer (optional)
AB = H H: Overload in checkweighing Y,Y,Y = Draft shield doors

A = L: Underweight XXX = Decimal value calculated from

AB = L L: Underweight in checkweighing binary data

Decimal value	Binary value	Control information
1	Bit 0 = 0:	No error/ionizer off
ı	Bit0 = 1:	Draft shield error/ionizer on
2	Bit $1 = 0$:	Draft shield motors off
Z	Bit $1 = 1$:	Draft shield in motion
0	Bit $3 = 0$:	Learning function off
8	Bit $3 = 1$:	Learning function on
16	Bit $4 = 0$:	At least one draft shield door open
10	Bit 4 = 1:	All draft shield doors closed
32	Bit $6 = 0$:	Motorized draft shield operation
ンム	D:4C 1.	Manual dueft shield an austion

R,M,L = COO: Right door closed(Closed), Middle and Left doors open (Open) R,M,L = OCC: Right door open (Open), Mmiddle and Left doors closed (Closed)

Error Message

Position 1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16
*	*	*	Е	r	r	*	*/#	#	#	*	*	*	*	CR	LF

^{*:} Spaces # # #: Error code number

Example: Output of the weight value + 111.255 g

Bit 6 = 1:

Position	1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	
	*	*	*	1	1	1		2	5	5	*	g	*	*	CR	LF	

Position 1: Plus +, or minus – or space Position 2: Space or weight value

Position 3 – 10: Weight value with decimal point, leading zeros are output as

spaces.

Position 11: Spaces

Position 12 - 14: Characters for unit of measure or space

Position 15: Carriage return
Position 16: Line feed

Data Output Format with 22 Characters

When data is output in this format, ID codes with 6 characters will precede data with a 16-character format. These six characters identify the subsequent value.

Normal Operation

K: ID code character¹)

*: Spaces

A: Displayed characters

LF: Line feed

E: Unit symbol character¹)

See chapter "Mass Unit Conversion"

CR: Carriage return

Special Outputs

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18	19	20	21	22
S	t	a	t	*	*	*	*	*	*	*	*	-	-	*	*	*	*	*	*	CR	LF
												Н	Н								
												L	L								
												С									

*: Spaces

- -: Final readout

H: Overload

H H: Overload in checkweighing

L: Underweight

LL: Underweight in checkweighing

C: Adjustment

Draft shield and ionizer status similar to data output format with

16 characters

Error Message

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18	19	20	21	22
S	t	a	t	*	*	*	*	*	Е	r	r	*	#	#	#	*	*	*	*	CR	LF

^{*:} Spaces # # #: Error code number

ID Code the Character K



The ID code depends on the model, e. g. not all units and IDs are available for verified balances

Device Information, Basic Data

Manuf.	Manufacturer
Model	Model
SerNo.	Serial number
BAC:	Balance version
APC:	Display version

Print function

Name Name

Task profiles

Task	Task title
TskDe	s Task description

User profiles

User User title
UsrDes User description

Data Memory

MemNo. Memory number
MemID Memory ID

Basic Weighing

G#	Gross weight 2	
N	Net weight	
N 1	Net weight (T1·0)	
T	Tare weight	
T1	Tare weight 1	
T 2	Tare weight 2	

Calibration/Adjustment

Target weight
Status
Deviation weight
Result ReproTEST

Leveling

LevRes Leveling result

SQmin Function

SQmin Minimum weight

Individual Identifiers

P-ID1	Print ID1
T-ID1	Task start ID1
I-ID1	Initialization ID1
R-ID1	Result output ID1
C-ID1	Component ID1
E-ID1	Evaluation ID1
X-ID1	Task end ID1

Density Determination

Density De	cermination
ID-Dk	Name of density determination kit
n	No. of wires
d	Wire diameter
D	Vessel diameter
Id-Liq	Name of liquid
RefFl	Density of liquid
b	Coefficient of expansion of the liquid
LA	Correction for air density
Pl.vol	Volume of glass plummet
Temp	Temperature of liquid
RhoFl	Density of liquid
Corr	Correction factor of the density determination kit
Wa	Weight of sample in air
Wfl	Weight of sample in liquid or weight of liquid
Wt	Residual liquid
Wr	Sample plus weight of liquid
Rho	Density of sample
Vol	Volume of sample

Statistics

n D e f	Default number of items
n	Transaction counter
x	Average value of the weight values
s	Standard deviation of the weight values
sRel	Variation coefficient of the weight values
Sum	Total of all weight values
Max	Highest weight value
Min	Lowest weight value
Dif	Difference between highest and lowest weight value

ID Code the Character K
(Continued)

Calculation

Form	Formula
Res	Results using equations
Χ	Variable X
Υ	Variable Y
а	Variable a
b	Variable b

Animal Weighing

mDef	Target value
Mul	Calculation factor
mAct	Actual number
x-Net	Results
x-Res	Calculated result

Formulation

i oi iiiaiati	UII	
IDCm1	ld Component 1	
TCmp1	Reference weight Comp. 1	
PCmp1	Reference percentage weight	
	Comp. 1	
Comp1	Weight value Comp.1	
WD1	Difference Weight 1	
PD1	Difference Percent 1	
WghMod Mode for inputting initial		
	weight	
n D e f	Default number	
	of components	
Total	Total of weight values	
RSum	Total reference	
n	Transaction counter	
IniWgt	Default total weight	

Weighing in Percent

pRef	Percentage of reference
Wxx%	Reference percentage weight
Prc	Percent
D	Percent loss
DR	Ratio in percent 1
0 R	Ratio in percent 2

Timer-controlled Functions

Tillici -coi	iti olicu i ulictiolis
InvTim Interval	
FixTim Fixed point in time	
ResTim Remaining time	
Hold	Display blocked

Totalizing

n D e f	Default number of items
Comp	Weight value start
Comp	Current weight value
n	Number of the item
CompC	Weight value start:
	calculated components
CompC	Current weight value:
	calculated components
Sum	Total of all weight values

DKD Measurement Uncertainty

a 1	Addend a1
b1	Factor b1
b1 e-	Exponent b1
a 2	Addend a2
b 2	Factor b2
b2 e-	Exponent b2
P fact	Pa Factor
P fact U	Pa Factor absolute Inaccuracy of the
	absolute Inaccuracy of the
U	absolute Inaccuracy of the process
U	absolute Inaccuracy of the process relative Inaccuracy of the

Parts Counting

wRef	Reference piece weight
n R e f	Reference sample quantity
Qnt	Piece count

Checkweighing

Setp	Target value
Min	Min. tolerance for checkw.
Max	Upper tolerance for checkw.
Lim-	Lower percentage
Lim+	Upper percentage
Count	Counter for correct values
CountLL	Counter for too small values
CountHH	Counter for too large values
Diff	Check result measure
	·-

Motorized Draft Shield

WPC	Draft shield version
Stat	Draft shield status
lonizer	

. .

Stat lonizer status

Data Input

SBI Commands (Data Input Format)

The computer connected via the data port can send commands to the balance to control balance and application program functions.

These control commands may have different formats and contain up to 26 characters. Each of these characters must be sent based on the setup configuration for data transmission.

Formats for Control Commands

Format 1:	Esc	!	CR	LF	
Format 2:	Esc	!	#	_	CR LF
Format 3:	Esc	!	#	Et	(max. 20 &) & _ CR LF

Esc: Escape

!: Command character

#: Number

8t: Number or letter_: Underline (ASCII: 95)CR: Carriage return (optional)LF: Line feed (optional)

max: Depends on command character (parameters): The entry is truncated after the max. length, and not rejected as when entered via the keyboard.

Examples:

Format 1: Esc P
Format 2: Esc x1_
Format 3: Esc t Article1_
Format 4: Esc e!{MESSAGE.SHOW_ERR

Format 4: Esc e!{MESSAGE.SHOW_ERROR "Weight too low!}_')
Esc e?{WEIGH.WGT_NET}_')

(e.g. query net weight value)

Esc ea{Art.413}_1)

(determine weight and tare values and save to the alibi memory together with the label)

1) Additional parameters and module identifications are available in the document "SBI Interface" as a download from the Sartorius homepage.

Overview of SBI Commands

Format	Command	Action/Function		Comments
1	ESC P	Print to the interface where the pro	mpt originated	Corresp. to Menu, with/without stability
1	ESC T	"TARE" key taring and zeroing		
1	ESC K	Filter "Very stable conditions"		
1	ESC L	Filter "Stable conditions"		
1	ESC M	Filter "Unstable conditions"		
1	ESC N	Filter "Very unstable conditions"		
1	ESC 0	Block keys		
1	ESC Q	Acoustic signal		
1	ESC R	Unblock keys		
1	ESC S	Restart		
1	ESC Z	Internal calibration/adjustment		Depending on menu, 1/2 step increments
1	ESC U	Tare		
1	ESC V	Zero key		
1	ESC W	Ext. Adjustment with default weight	t	Depending on menu, 1/2 step increments
2	ESC f3_	Zero		
2	ESC f4_	Tare (without zeroing)		
2	ESC f5_	Left draft shield key (closing and op	ening as learned or default)	Only if available
2	ESC f6_	Right draft shield key(closing and or	pening as learned or default)	Only if available
2	ESC kF1_	Soft key 1, first from right		Depends on application
2	ESC kF2_	Soft key 2, second from right		Depends on application
2	ESC kF3_	Soft key 3, third from right		Depends on application
2	ESC kF4_	Soft key 4, fourth from right		Depends on application
2	ESC kF5_	Soft key 5, fifth from right		Depends on application
2	ESC kF7_	"Menu" soft key		
2	ESC kF9_	"TASK" key		
2	ESC kF10_	"USER" key		
2	ESC kP_	Print as with "PRINT" key (e. g. to so	everal interfaces)	
2	ESC m0_	lonizer status		Only if available
2	ESC m1_	lonizer on, with preset time		Only if available
2	ESC m2_	lonizer off		Only if available
2	ESC s3_	Back, exit, cancel		
		Models with the	Models with the	Analytical and
	TCO o	analytical draft shield	rotational draft shield	rotational draft shield
2	ESC w0_	Draft shield status	Draft shield status	Only if available
2	ESC w1_	Open left door Close all doors	Open draft shield 100% to the left Close draft shield	Only if available
2	ESC w2_ ESC w3_			Only if available Only if available
2		Open upper door	Open draft shield door 100% to the right	<u> </u>
$\frac{2}{2}$	ESC w4_ ESC w5_	Open right door	Open draft shield door 100% to the right	Only if available Only if available
2		Open left and upper doors Open left and right doors	_	Only if available
2	ESC w6_ ESC w7_	Open right and upper door		Only if available
$\frac{2}{2}$	ESC w8_	Open all doors	-	Only if available
2	ESC x0_	Perform internal calibration		Offig it available
2	ESC x1_	Print weigher type		
2	ESC x2	Print serial number		
2	ESC x3_	Print balance software version		
2	ESC x4_	Print softw. version of display and c	ontrol unit	
2	ESC x5_	Print user/device 1D		
2	ESC x6_	Print cal./adjust. weight ID		
2	ESC x7_	Print batch ID		
3	ESC txxx_	Input text		
3		Enter opening angle xxx in degrees	for left key	
			•	
3		Enter opening angle xxx in degrees	tor ngnt key	
3	ESC z5xxx_	Input user/device ID		
3	ESC z6xxx_	Input cal./adjust. weight ID		
3	ESC z7xxx_	Input batch/lot ID)	
4	ESC e! {xxx}_	Execute application command, {xxx	,	
4	ESC e? {xxx}_	Query parameter, $\{xxx\}$ = Paramete		
4	ESC ea {xxx}_	Save weight value to alibi memory (v	with optional label), $\{xx\}$ = Memory ID	

SICS Commands

The computer connected via the data port can send control commands to the balance to control balance and application program functions.



Please consult the separate SICS documentation for a description of the syntax.

Command Action/Function

Level 0 Version 2.3x

- @ Reset all SICS commands
- I 0 List all available commands
- I 1 Send information about the level and its versions
- I 2 Balance model prompt
- I 3 Prompt for input of software version of balance (BAC)
- I 4 Prompt for balance serial number
- I 5 Prompt for software version of display (APC)
- **S** Send weight value at stability
- SI Send weight value without stability
- SIR Send automatic weight values at and without stability
 - Z Zero the balance at stability
 - Z I Zero the balance without stability

Level 1 Version 2.2x

- T Tare the balance at stability
- TI Tare the balance without stability
- **D** Write text in display
- **DW** Delete text from display
- K Key control
- TA Prompt for and allocate tare memory
- TAC Delete tare memory
- SR Send weight value if weight changes

Level 2

- M13 Activate/deactivate touch screen softkeys
- PWR Turn balance on/off (standby)
 - **SU** Send weight value at stability with current weight unit (with motorized draft shield)
- WS Query door position of motorized draft shield / Open or close door
- MO7 Query / Activate / Deactivate automatic draft shield
- M24 "Print" key query / deactivate / print the stable or unstable weight values

Remote control (remote)

- P112 Write text in selected line on display
- P113 Delete text from selected line in display
- P114 Overwrite task or user names
- P120 Turn off bar graph in checkweigher
- P121 Turn on bar graph in checkweigher
- RM20 Activate/deactivate user input
- **RM30** Assign new function to softkeys
- RM36 Assign/prompt for function of multiple softkey lines
- RM38 Activate RM36-assigned softkey lines
- RM39 Activate/deactivate RM30-assigned softkey functions

Additional Sartorius commands

- S A Send weight value at stability and store in alibi memory (with optional label)
- **CMD** Execute application command
- PAR Query parameter
- MN36 Assign a description to several menus
- MN38 Display or hide assigned menus assigned with MN36
- **TX36** Assign texts to several text pages
- TX37 Overwrite a line of a text page
- TX38 Activate / Deactivate assigned text pages with TX38

Updating the Software

Purpose:

To better serve our customers, Sartorius continues to improve its balance software. Sartorius makes the latest software versions available online so that you can quickly and easily remain up-to-date. Sartorius, however, makes no guarantee regarding the use of this software.

The software should only be updated by an administrator.




On verified balances, observe the Audit Trail instructions under (3) Uploading the Software Update to the Balance.



Before updating the software, save all data (see chapter Exporting Data).



If you are working with User Management (user profiles are defined), the software can only be updated by the administrator.

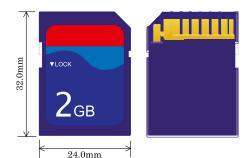


1) Download the software update from the Internet


The first step is to download the software to your computer:

- In your browser, go to website
 - "www.sartorius.com/downloads" and click on the "Software" link.
- Select the update packet for your balance.
- ► Enter the information required for registration.
- ▶ Download the software packet to your computer.
- Unpack the zip file.
- \triangleright Copy the software update to a SD card (32 × 24 × 2.1 mm).

2) Insert the memory card into the control unit

The slot for the SD card is located at the back of the control unit.

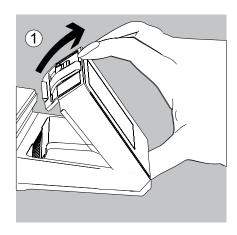

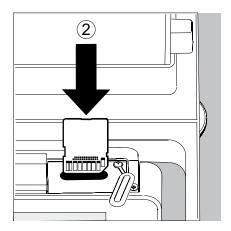

- 1. Tilt the display and control unit.
- 2. Swivel the cover of the card slot to the front.
 Align the SD card so that the contacts are facing forward.
- 3. Insert the card into the slot as far as it will go.
- ► To remove the card, press it against the resistance in the direction of the slot so that the card springs out.



3) Upload the software update to the balance:

- If you are working with User Management, it is necessary to log on as administrator.
- ► Toggle to the System Settings menu.
- ▶ Open the **Update software** submenu.
- > You are asked if you want to delete the existing profiles.
- ► Touch No or Yes.

Once the SD card with the current software is inserted, touch Next.

If the following error message appears:

Wrong XML version. Please check the data.

- ► Touch Next.
- ➤ All previous data will be saved in the new format.
- After updating your software, check your settings (Menu, TASK and USER).
- ▶ If the XML files continue to cause problems, contact your local Sartorius service center.
- ▶ After a version check is run, the versions of the current and the new software are displayed.
- To start the update, press Continue.
- An acoustical signal will sound when the update is complete and the balance then carries out a self test.

Audit Trail for Software Download on Models Verified for Legal Metrology

Software Status of the Balance

M

Continue

Models verified for use in legal metrology are assigned an authorized software status when the balance is shipped for delivery. This can be indicated in the display. If something concerning the software or the legal regulations changes, the software status must be updated.

Display Audit Trail of the Balance Software

- ► Turn off the balance and disconnect it from the power supply.
- Keep () pressed and reconnect the balance back to the power supply.
- ➤ The "Bios" menu is displayed.
- Press the "Settings" menu item.
- ▶ Select the "Audit trail of software updates" menu item.
- Select "Show audit trail."
- ➤ The Audit Trail with modification date, name, new and old software and version number are indicated in the display.

Back

Error and Status Messages

Error messages appear directly in the respective active application, usually with a descriptive text on how to correct it. Warning and reminder functions are displayed with a fixed 3-level action hierarchy:




Status display (tab):

The available status displays and their meanings are described in the "User interface" section.



Warning text with a repeating reminder

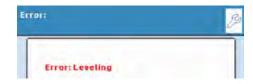

Error text with mandatory prompt for troubleshooting

You cannot proceed until you have eliminated the error.

Ineffective key commands or entries

If you press a key that has no function or not allowed, this error is indicated as follows:

- A double beep indicates that the key has no function assigned to it.
- An error message indicates invalid input.
- An error code or error message indicates incorrect operation.

Error messages are shown on the display for 2 seconds or must be acknowledged by pressing a key. The error text provides direct suggestions on how to proceed. If the error cannot be resolved by following the suggestions, please contact Sartorius customer service.

GPL License

Devices of the Cubis series also contain "free software" in the firmware that is licensed under the GNU General Public License (GPL) Version 2, June 1991 and the GNU Lesser General Public License (LGPL) Version 2.1, February 1999. This third-party developed "free software" is copyrighted and is provided free of charge.

The licensing conditions of the Free Software Foundation, Inc. are available in English on the included CD-ROM.

You can purchase the GPL source text from Sartorius on CD under the VF no. 4043 for a shipping cost of € 20.

Care and Maintenance

Service

Regular servicing by a Sartorius technician will extend the service life of your balance and ensure its continued weighing accuracy. Sartorius offers its customers service contracts with regular maintenance intervals ranging from 1 month to 2 years. The frequency of the maintenance intervals depends on the operating conditions and user's tolerance requirements.

Repairs

Repair work must only be performed by trained service technicians. Repairs performed by untrained persons may result in considerable hazards for the user.



The device should be unplugged during repair work.

Unplug the power cord from the outlet.

Repair work must only be performed by Sartorius-trained service technicians. As needed: Please contact your Sartorius dealer!


Cleaning the Balance




Ensure that no dust or liquid enters the balance weighing system.

Do not use aggressive cleaning agents (solvents or similar agents).


lsolate from supply voltage: Unplug the power cord from the wall outlet and unplug any data cables connected to the balance.

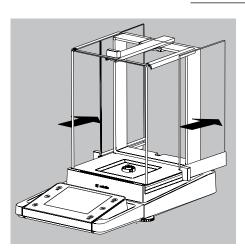


- ▶ Open the panels of the analytical draft shield completely.
- ► Clean the housing and interior of the balance with a cloth lightly moistened with a soap solution.
- Dry all parts with a soft, dry cloth or use blotting paper to absorb dampness. Then replace all dried parts.
- ▶ Remove panels from the analytical draft shield and clean them with commercially available glass cleaner. Dry the panels and then reinstall them in the balance.

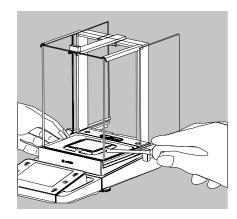

Carefully remove any sample residue/spilled powder using a brush or hand-held vacuum.

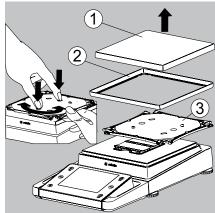

- ▶ If necessary, remove the weighing pan, shield plate and pan support.
- 1) Weighing pan
- 2) Shield plate/draft shield
- 3) Pan support

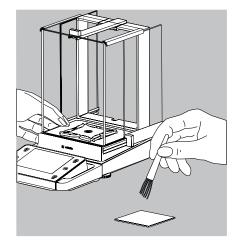

► Clean parts with a cloth or brush.

Then replace the parts.

Cleaning Stainless Steel Surfaces

All stainless steel parts should be cleaned at regular intervals.

Use a damp cloth or sponge to clean stainless steel parts on the balance. Only use conventional household cleaning agents which are suitable for stainless steel (e.g. Stahlfix). After this, let the device dry.


For additional protection, protective oil may be applied.

Remove the stainless steel weighing pan and thoroughly clean it separately. Only use solvents for cleaning stainless steel parts.

The stainless steel weighing pan should be cleaned simply by rubbing. After this, rinse the equipment thoroughly until all residue is removed.

No protective oil should be applied to the stainless steel weighing pan!

Safety Inspection

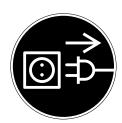

If there is any indication that safe operation of the balance is no longer warranted:

- Disconnect from the supply voltage: Unplug the power cord from the outlet.
- Secure the AC adapter and cord so that they cannot be used.

Safe operation of the mass comparator with the AC adapter is no longer ensured when:

- The AC adapter or the mains connecting lead shows visible damage.
- The AC adapter no longer functions properly.
- Following extended storage in adverse conditions.
 In this case, notify the Sartorius Service Center.

Maintenance and repair work may be performed only by authorized service technicians who have access to the required maintenance manuals and instructions and who have received the necessary training.

We recommend that the AC adapter be inspected by a qualified service technician with regard to the following:

- Leakage current: <0.25 mA measured with a properly calibrated multimeter.
- Insulation resistance >7M0hm as measured with a constant voltage of at least 500 volts at a 500 k0hm load.


The duration and number of checks should be determined by a qualified Sartorius service technician on site based on specific ambient and operating conditions (once a year as a minimum).

Shipping the Balance


If repairs are required, use the original packaging to transport the balance. To ensure adequate protection for safe shipment, Sartorius products have been packaged to the extent necessary using environmentally friendly materials. Only the original packaging provides optimum protection for the equipment.




- Disconnect the device from the power supply.
- Disconnect any data cables from the device.

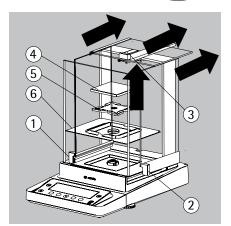


- ▶ Remove all items (such as weights, sensors, etc.) from the weighing chamber.
- 1) Remove the other side panel
- **2)** Remove the other side panel
- **3)** Remove the upper draft shield panel
- **4)** Remove the weighing pan
- **5)** Remove the pan support
- 6) Remove the shield plate/draft shield

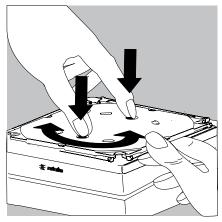

- ▶ On models without a draft shield: press down on the two pan support fasteners.
- ▶ Rotate and remove the pan support.

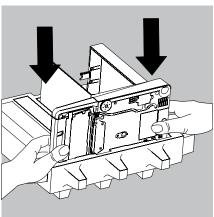


▶ Place the balance in the lower part of the packaging.

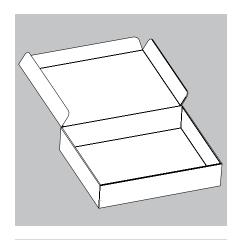


Transporting the Parts (Large Analytical Draft Shield)


▶ Get the box for the individual parts of the balance ready.

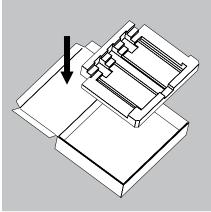


▶ Place the bottom foam piece in the box.

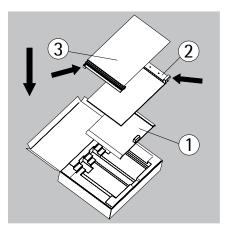

- ▶ Place the panels in the packaging:
- 1) Place the upper draft shield panel into the packaging (handle upwards).
- 2) Place the side draft shield panel into the packaging (handle upwards).
- **3)** Place the other side panel into the packaging (handle downwards).

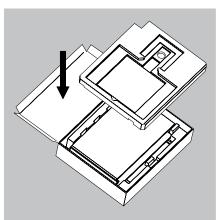


► Then place the top foam piece in the box.

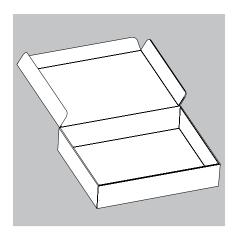


Only for weighing pan:

Place the shield plate into the box.

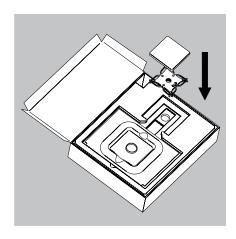


- ▶ Place the pan support and weighing pan into the opening.
- Close the box.

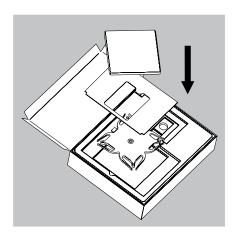


Only for large weighing pan:

- ▶ Place the following parts into the foam piece in the order given:
- 1) Pan support
- 2) Shield plate
- 3) Weighing pan
- Close the box.

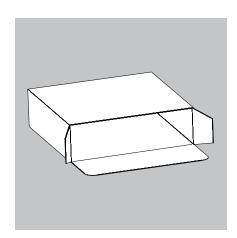


Transporting the Parts (Small Analytical Draft Shield)


► Get the box ready for the individual parts of the balance.

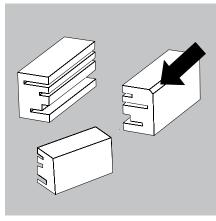


► Get the foam pieces ready.

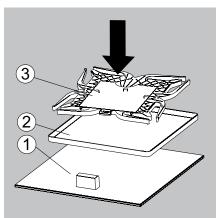


- Place the balance parts on top of each other.
- 1) Upper draft shield panel
- 2) Shield plate/draft shield
- 3) Pan support

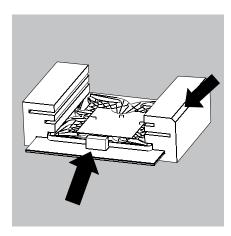


Slide the parts into the foam.

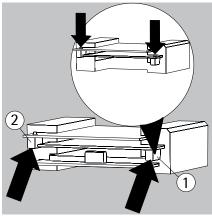


- ▶ Place the panels in the packaging.
- 1) Place the side panel into the packaging (handle downwards).
- 2) Place the side panel into the packaging (handle downwards).

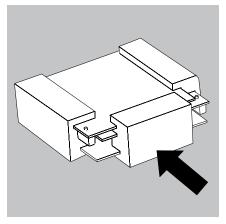


▶ Place the foam piece in front of the parts.

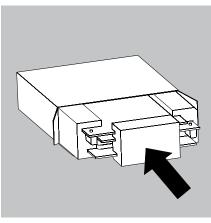


▶ Place the package into the box.

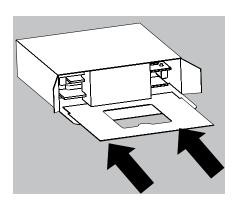


Slide the shield plate into the packaging.

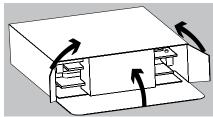


Close the box.

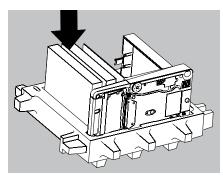

Place the box into the packaging.

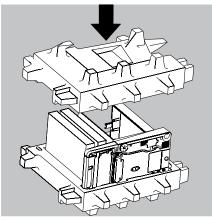

Place the top part on to the packaging.

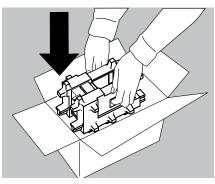

Insert the balance into the box with cushioning.

Send the packaged balance.

Disposal



The packaging is to be taken to a local waste disposal site if no longer required. The packaging comprises of environmentally-friendly materials that can be used as secondary raw materials. The equipment, including accessories and batteries, should not be disposed of as regular household waste. EU legislation in Member States requires electrical and electronic equipment to be collected separately from unsorted municipal waste so that it may be recycled. In Germany and several other countries, Sartorius itself assumes responsibility for the return and conformant disposal of its electronic and electrical products. These products may not be placed with household waste or brought to collection centers run by local public disposal operations – not even by small commercial operators. For disposal in Germany and in the other member nations of the European Economic Area, please contact our local service technicians or our Service Center in Goettingen:

Sartorius Service Center, Weender Landstrasse 94-108, 37075 Goettingen, Germany

In countries that are not members of the European Economic Area (EEA) or where no Sartorius subsidiaries or dealerships are located, please contact your local authorities or a commercial disposal operator. Prior to disposal and |or scrapping of the equipment, any batteries should be removed and disposed of in local collection boxes. Sartorius will not take back equipment contaminated with hazardous materials (ABC contamination) – either for repair or disposal.

Please refer to our web site (www.sartorius.com) or contact the Sartorius Service Department for more detailed information regarding repair service addresses or the disposal of your device.

Specifications

General Data

Primary	100 – 240V~, -15%/+10%, 50-60Hz, 1.0A
Secondary	15V, \pm 5%, 2.66A (max.), protected electronically against short circuit
Power supply connection cable	Two-sided plug with a 3-pin country-specific power plug and 3-pin socket (IEC/EN60320-1/C14) for connection to the power supply
Other data	See label on the power supply or the user manual on
Balance	
Power supply	Only via Sartorius power supply 6971987
Input voltage	15 VDC, ± 5%
Power consumption	15 W (max.)
Ambient conditions	
Environment	For indoor use only
Ambient temperature: Storage and shipping	−10 °C to +60 °C
Ambient temperature: Operation	+5 °C to +40 °C
Height	2000 m above sea level
Highest relative humidity: humidity for 40 °C	80% for temperatures up to 31 °C, decreasing linearly up to 50% relative
Safety of electrical equipment	According to EN 61010-1:2001 Safety requirements for electrical equipment for measurement, control, and laboratory use - Part 1: General requirements
Electromagnetic compatibility Defined immunity to interference: Limitation of Emissions:	In accordance with EN 61326-1:2006 Electrical equipment for measurement, control, and laboratory use EMC requirements – Part 1: General requirements Suitable for use in industrial areas Class B (suitable for use in residential areas and areas that are connected
Standard equipment	to a low voltage network that also supplies residential buildings).
Customization to operating	
and setup conditions	By selection of 1 of 4 optimized filter levels
Selectable weight units	Gram, Kilogram, Carat, Pound, Ounce, Troy Ounce, Tael Hong Kong, Tael Singapore, Tael Taiwan, Grain, Pennyweights, Milligram, Parts per Pound, Tael China, Mommes, Austrian carat, Tola, Baht and Mesgh
Available application programs	Changing unit, counting, weighing in percent, animal weighing, calculation, density determination, checkweighing, timer controlled functions, totalizing, formulation, statistics, 2nd tare memory, identifiers, product data memory

Microbalances 0.001mg

Model		MSA6.6S	MSA6.6S-F	MSA3.6P
Readability	mg	0.001	0.001	0.001/0.002/0.005
Weighing capacity	g	6.1	6.1	1.1/2.1/3.1
Tare range (subtractive)	g	-6.1	-6.1	-3.1
Repeatability	≤±mg	0.001	0.001	0.003/0.004/0.005
Linearity	≤±mg	0.004	0.004	0.004
Corner load (test load [g])	mg	4 (2 g)	4 (2 g)*	5 (1 g)
Min. initial weight	mg	2	-	4
Sensitivity drift between +10+30°C	±ppm/K	1	1	1
Typical stabilization time	S	≤ 5	≤ 5	≤ 5
Typical measurement time	S	≤ 8	≤ 8	≤ 8
External standard calibration value (min. accuracy class)	g	5 (E2)	5 (E2)	3 (E2)
Display result				
(depending on the set filter level)		0.1 - 0.4	0.1 - 0.4	0.1 - 0.4
Weighing pan size \varnothing	mm	30	50	30
Weighing chamber height	mm	70	-	70
Protection		Protected against dust	and water	

Ultra-Microbalances 0.0001mg

Model		MSA2.7S	MSA2.7S-F
Readability	mg	0,0001	0,0001
Weighing capacity	g	2.1	2.1
Tare range (subtractive)	g	-2.1	-2.1
Repeatability	≤±mg	0.00025	0.00025
Linearity	≤±mg	0.0009	0.0009
Corner load (test load [g])1)	mg	2.5 (1)	2.5 (1)
Min. initial weight	mg	1	-
Sensitivity drift between zw. +10+30°C	±ppm/K	1	1
Typical stabilization time	S	≤ 7	
Typical measurement time	S	≤ 10	
External standard calibration value (min. accuracy class)	g	2 (E2)	2 (E2)
Display result			
(depending on the set filter level)		0.1 - 0.4	0.1 - 0.4
Weighing pan size \varnothing	mm	20	50
Weighing chamber height	mm	70	15
Protection		Protected against dust a	and water

¹) = Standard pan

Semi-microbalances 0.01 mg

Model		MSA225S	MSA225P	MSA125P
Readability	mg	0.01	0.01/0.02/0.05	0.01/0.1
Weighing capacity	g	220	60/120/220	60/120
Tare range (subtractive)	g	- 220	- 220	- 120
Repeatability	≤±mg	0 to 6 g: 0.015 60 to 220g: 0.025	0 to 6 g: 0.015 60 to 220g: 0.04	0 to 6 g: 0.015 60 to 120g: 0.06
Linearity	≤±mg	0.1	0.15	0.15
Corner load (test load [g])	mg	0.15 (100)	0.2 (100)	0.15 (50)
Min. initial weight*	mg	20	20	20
Sensitivity drift between +10 to +30°C	±ppm/l	K 1	1	1
Typical stabilization time	S	≤ 2	≤ 2	≤ 2
Typical measurement time	S	≤ 6	≤ 6	≤ 6
External standard calibration value (min. accuracy class)	g	200 (E2)	200 (E2)	100 (E2)
Display result (depending on the set filter level)		0.2 - 0.4		
Weighing pan size (W × D)	mm	85 x 85		
Weighing chamber height (draft shield DU)	mm	261		
Protection		Protected against dust	and water	

^{* =} Typical min. initial weighing according to USP (Unites States Pharmacopeia), USP31-NF26

Analytical balances 0.1 mg

Model		MSA524S	MSA524P	MSA324S	MSA224S	MSA324P	MSA124S
Readability	mg	0.1	0.1/0.2/0.5	0.1	0.1	0.1/0.2/0.5	0.1
Weighing capacity	g	520	120/240/520	320	220	80/160/320	120
Tare range (subtractive)	g	- 520	- 520	- 320	- 220	- 320	- 120
Repeatability	<±mg	0.1	0.15/0.2/0.4	0.1	0.07	0.1/0.2/0.4	0.1
Linearity	<±mg	0.4	0.5	0.3	0.2	0.5	0.2
Corner load (test load [g])	mg	0.3 (200)	0.4 (200)	0.3 (200)	0.2 (100)	0.4 (200)	0.2 (50)
Min. initial weight*	mg			120	120	120	120
Sensitivity drift between +10 to +30°C	±ppm/	K 1	1	1	1	1	1
Typical stabilization time	S	< 1	< 1	< 1	< 1	< 1	< 1
Typical measurement time	S	< 3	< 3	< 3	< 3	< 3	< 3
External standard calibration value (min. accuracy class)	g	500	500	200+100 (E2)	200 (E2)	200+100 (E2)	100 (E2)
Display result (depending on the set filter level)		0.1 - 0.4					
Weighing pan size (W × D)	mm	85 × 85					
Weighing chamber height (draft shield DU)	mm	261					
Protection		Protected a	gainst dust an	ıd water			

^{* =} Position according to OIML R76

Precision balances

Models		MSA5203S	MSA5203P	MSA3203S	MSA3203P		
Readability	mg	1	1/2/5	1	1/10		
Weighing capacity	g	5200	1200/2400/5200	3200	1010/3200		
Tare range (subtractive)	g	- 5200	- 5200	- 3200	- 3200		
Repeatability	≤±mg	1	1	1	1/6		
Linearity	≤±mg	5	5	5	5		
Corner load (test load [g])	mg	2 (2000)	2 (2000)	2 (1000)	2 (1000)		
Min. initial weight*	g				1.5		
Sensitivity drift between +10 to +30°C	±ppm/l	K 1	1	1	1		
Typical stabilization time	S	≤ 1	≤ 1	≤ 1	≤ 1		
Typical measurement time	S	≤ 2	≤ 2	≤ 2	≤ 1.5		
External standard calibration value (min. accuracy class)	g	5000	5000	2000	2000 (E2)		
Display result (depending on the set filter level)		0.1 - 0.4					
Weighing pan size (W × D)	mm	140 × 140					
Weighing chamber height (draft shield DE)	mm	172					
Protection		Protected against dust and water					

^{* =} Position according to OIML R76

Models		MSA2203S	MSA2203P	MSA1203S
Readability	mg	1	1/10	1
Weighing capacity	g	2200	1010/2200	1200
Tare range (subtractive)	g	- 2200	- 2200	- 1200
Repeatability	≤±mg	1	1/6	0.7
Linearity	≤±mg	3	5	2
Corner load (test load [g])	mg	2 (1000)	3 (1000)	2 (500)
Min. initial weight*	g	1.5	1.5	1.5
Sensitivity drift between +10 to +30°C	±ppm/	K 1	1	1.5
Typical stabilization time	S	≤ 1	≤ 1	≤ 1
Typical measurement time	S	≤ 1.5	≤ 1.5	≤ 1.5
External standard calibration value (min. accuracy class)	g	2000 (E2)	1000 (E2)	1000 (E2)
Display result (depending on the set filter level)		0.1 - 0.4		
Weighing pan size (W \times D)	mm	140 × 140		
Weighing chamber height				
(draft shield DE)	mm	172		
Protection		Protected agains	t dust and water	

^{* =} Position according to OIML R76

Precision balances

Models		MSA623S	MSA623P	MSA323S
Readability	mg	1	1/2/5	1
Weighing capacity	g	620	150/300/620	320
Tare range (subtractive)	g	- 620	- 620	- 320
Repeatability	≤±mg	0.7	1/2/4	0.7
Linearity	≤±mg	2	5	2
Corner load (test load [g])	mg	2 (200)	4 (200)	2 (200)
Min. initial weight*	g	1.5	1.5	1.5
Sensitivity drift between +10 to +30°C	±ppm/	K2	2	2
Typical stabilization time	S	≤ 0.8	≤ 0.8	≤ 0.8
Typical measurement time	S	≤ 1	≤ 1	≤ 1
External standard calibration value (min. accuracy class)	g	500 (E2)	500 (F1)	200 (E2)
Display result (depending on the set filter level)		0.1 - 0.4		
Weighing pan size (W × D)	mm	140 × 140		
Weighing chamber height (draft shield DE)	mm	172		
Protection		Protected agains	t dust and water	

^{* =} Position according to OIML R76

Models		MSA14202S	MSA14202P	MSA10202S	MSA8202S	
Readability	mg	10	10/20/50	10	10	
Weighing capacity	g	14,200	3500/7000/14,200	10,200	8200	
Tare range (subtractive)	g	- 14,200	- 14,200	- 10,200	- 8200	
Repeatability	≤±mg	10	10/20/40	7	7	
Linearity	≤±mg	30	50	20	20	
Corner load (test load [g])	mg	20 (5000)	40 (5000)	20 (5000)	20 (5000)	
Min. initial weight*	g	15	15	12	12	
Sensitivity drift between +10 to +30°C	±ppm/	K 1.5	1.5	2	2	
Typical stabilization time	S	≤ 1	≤ 1	≤ 1	≤ 1	
Typical measurement time	S	≤ 1.5	≤ 1.5	≤ 1.5	≤ 1.5	
External standard calibration value (min. accuracy class)	kg	10 (E2)	10 (E2)	10 (E2)	5 (E2)	
Display result						
(depending on the set filter level)		0.1 - 0.4				
Weighing pan size $(W \times D)$	mm	206×206				
Protection		Protected against dust and water				

^{* =} Position according to OIML R76

Precision balances

Models		MSA6202S	MSA6202P	MSA5202S	MSA4202S
Readability	mg	10	10/20/50	10	10
Weighing capacity	g	6200	1500/3000/ 6200	5200	4200
Tare range (subtractive)	g	- 6200	- 6200	- 5200	- 4200
Repeatability	≤±mg	7	7/20/40	6	7
Linearity	≤±mg	20	50	10	20
Corner load (test load [g])	mg	20 (2000)	50 (2000)	10 (2000)	30 (2000)
Min. initial weight*	g	12	12	10	12
Sensitivity drift between +10 to +30°C	±ppm/	K 2	2	2	2
Typical stabilization time	S	≤ 1	≤ 1	≤ 0.8	≤ 0.8
Typical measurement time	S	≤≤ 1.5	≤ 1.5	≤ 1	≤ 1
External standard calibration value (min. accuracy class)	kg	5 (E2)	5 (F1)	5	2 (E2)
Display result					
(depending on the set filter level)		0.1 - 0.4			
Weighing pan size $(W \times D)$	mm	206 × 206	206×206	140 × 140	206×206
Protection	·	Protected again	nst dust and water		

^{* =} Position according to OIML R76

Models		MSA2202S	MSA1202S	MSA12201S	MSA8201S	MSA5201S
Readability	mg	10	10	100	100	100
Weighing capacity	g	2200	1200	12,200	8200	5200
Tare range (subtractive)	g	- 2200	- 1200	- 12,200	- 8200	- 5200
Repeatability	≤±mg	7	7	50	50	50
Linearity	≤±mg	20	20	100	100	100
Corner load (test load [g])	mg	20 (1000)	20 (500)	200 (5000)	200 (5000)	200 (2000)
Min. initial weight*	g	12	12	100	100	100
Sensitivity drift between +10 to +30°C	±ppm/	K 2	2	4	4	4
Typical stabilization time	S	≤ 0.8	≤ 0.8	≤ 0.8	≤ 0.8	≤ 0.8
Typical measurement time	S	≤ 1	≤ 1	≤ 1	≤ 1	≤ 1
External standard calibration value (min. accuracy class)	kg	2 (F1)	1 (F1)	10 (F1)	5 (F2)	5 (E2)
Display result						
(depending on the set filter level)		0.1 - 0.4				
Weighing pan size $(W \times D)$	mm	206×206				
Protection		Protected again	nst dust and w	ater		

^{* =} Position according to OIML R76

Verified Models with EC Type Approval Certificate: Micro- and ultramicrobalances

Model		MSE6.6S-0CE	MSE2.7S-0CE	MSE3.6P-0CE	
Accuracy class*		I	I		
For verified models: EC Type Approval	Certificate	D09-09-015, Type	: MSX		
Scale interval d*	mg	0.001	0.0001	0.001	
Weighing capacity max*	g	6.1	2.1	3.1	
Calibration value e*	mg	1	1	1	
Min. load min*	mg	0.1	0.01	0.1	
Tare equalization range (subtractive)		≤ 100% from max	≤ 100% from max. weighing capacity		
Application range according to DIR*	g	0.001 - 6.1	0.001 - 2.1	0.001 - 3.1	
Min. initial weight**	mg	2	1	4	
Typical stabilization time	S	≤ 5	≤ 7	≤ 5	
Typical measurement time	S	≤ 8	≤ 10	≤ 8	
External standard calibration value (min. accuracy class)	g	5	2	3	
Application range (temperature)			nction: +5+40°C function: +15 +2	25°C	
Display result (depending on the set fil	ter level)	By selection of 1 of	of 4 optimized filter	levels	
Weighing pan size \varnothing	mm	30	20	30	
Weighing chamber height (draft shield DM)	mm	70	70	70	
Protection		Protected against	dust and water		

^{*} DIR = Directive 90/384/EEC on non-automatic weighing instruments used within the European Economic Area

** = Typical min. initial weighing according to USP (Unites States Pharmacopeia), USP31-NF26

Verified models with EC Type Approval Certificate: Semi-microbalances 0.01 mg

Model		MSA225S-0CE	MSA225P-0CE	MSA125P-0CE		
Accuracy class*		I	I			
For verified models: EC Type-Approval	Certificat	e D09-09-015, Type: 1	MSX			
Scale interval d*	mg	0.01	0.01/0.02/0.05	0.01/0.1		
Weighing capacity max*	g	220	60/120/220	60/120		
Calibration value e*	mg	1	1	1		
Min. load min*	mg	1	1	1		
Tare equalization range (subtractive)	≤ 100% from max. v	≤ 100% from max. weighing capacity				
Application range according to DIR*	g	0.001 - 220	0.001 - 220	0.001 - 120		
Min. initial weight**	mg	20	20	20		
Typical stabilization time	S	≤ 2	≤ 2	≤ 2		
Typical measurement time	S	≤ 6	≤ 6	≤ 6		
External standard calibration value (min. accuracy class)	g	200 (E2)	200 (E2)	100 (E2)		
Application range (temperature)		With "isoCAL" function	on: +5 to +40 °C Without "	isoCAL" function: +15 to +25 °C		
Adaptation to ambient conditions		By selection of 1 of 4 optimized filter levels				
Display result (depending on the set file	ter level)	0.2 - 0.4				
Weighing pan size (W × D)	mm	85 × 85				
Weighing chamber height (draft shield Di	J) mm	261				
Protection		Protected against du	ıst and water			

Verified models with EC Type Approval Certificate: Analytical balances 0.1 mg

Model		MSA524S- OCE	MSA524P- OCE	MSA324S- 0CE	MSA224S- OCE	MSA324P- 0CE	MSA124S- OCE
Accuracy class*		I	I	I	I	I	I
For verified models: EC Type-Approval Co	ertificate	D09-09-015	, Type: MSX				
Scale interval d*	mg	0.1	0.1/0.2/0.5	0.1	0.1	0.1/0.2/0.5	0,1
Weighing capacity max*	g	520	120/240/520	320	220	80/160/320	120
Calibration value e*	mg	1	1	1	1	1	1
Min. load min*	mg	10	10	10	10	10	10
Tare equalization range (subtractive)		≤ 100% from	n max. weighir	ng capacity			
Application range according to DIR*	g	0.01-520	0.01-520	0.01-320	0.01-220	0.01-320	0.01-120
Min. initial weight**	mg	120	120	120	120	120	120
Typical stabilization time	S	≤ 1	≤ 1	≤ 1	≤ 1	≤ 1	≤ 1
Typical measurement time	S	≤ 3	≤ 3	≤ 3	≤ 3	≤ 3	≤ 3
External standard calibration value (min. accuracy class)	g	500	500 (E2)	200+100 (E2)	200 (E2)	200 +100 (E2)	100
Application range (temperature)		With "isoCAL	" function: +5	to +40 °C Wi	thout "isoCAl	_" function: +1	5 to +25 °C
Display result (depending on the set filter	level)	0.1 - 0.4					
Weighing pan size (W × D)	mm	85 × 85					
Weighing chamber height (draft shield DU)	mm	261					
Protection		Protected ac	gainst dust and	d water			

^{*} DIR = Directive 90/384/EEC on non-automatic weighing instruments used within the European Economic Area

** = Typical min. initial weighing according to USP (Unites States Pharmacopeia), USP31-NF26

Verified models with EC Typ	Approval Certificate:	Precision balances
-----------------------------	-----------------------	--------------------

Models		MSA5203S-0CE	MSA5203P-0CE	MSA3203S-0CE	MSA3203P-0CE	
Accuracy class*		I		I	I	
For verified models: EC Type-Approval C	ertificate	D09-09-015, Type	: MSX			
Scale interval d*	mg	1	1/2/5	1	1/10	
Weighing capacity max*	g	5200	1200/2400/5200	3200	1010/3200	
Calibration value e*	mg	10	10	10	10	
Min. load min*	mg	100	100	100	100	
Tare equalization range (subtractive)		≤ 100% from max.	weighing capacity			
Application range according to DIR*	g	0.1 - 5200	0.1 - 5200	0.1 - 3200	0.1 - 3200	
Min. initial weight**	g	1.5	1.5	1.5	1.5	
Typical stabilization time	S	≤ 1	≤ 1	≤ 1	≤ 1	
Typical measurement time	S	≤ 2	≤ 2	≤ 2	≤ 1.5	
External standard calibration value (min. accuracy class)	g	5000	5000	2000	2000 (E2)	
Application range (temperature)		With "isoCAL" function: +5 +40 °C Without "isoCAL" function: +15 +25 °C				
Display result						
(depending on the set filter level)		0.1 - 0.4				
Weighing pan size (W \times D)	mm	140 × 140				
Weighing chamber height (draft shield DE)	mm	172				
Protection		Protected against of	dust and water			

Models		MSA2203S-0CE	MSA2203P-0CE	MSA1203S-0CE	
Accuracy class*		I	I		
For verified models: EC Type-Approval Co	ertificate	D09-09-015, Type:	MSX		
Scale interval d*	mg	1	1/10	1	
Weighing capacity max*	g	2200	1010/2200	1200	
Calibration value e*	mg	10	10	10	
Min. load min*	mg	100	100	100	
Tare equalization range (subtractive)		≤ 100% from max. v	weighing capacity		
Application range according to DIR*	g	0.1 - 2200	0.1 - 2200	0.1 - 1200	
Min. initial weight**	g	1.5	1.5	1.5	
Typical stabilization time	S	≤ 1	≤ 1	≤ 1	
Typical measurement time	S	≤ 1.5	≤ 1.5	≤ 1.5	
External standard calibration value (min. accuracy class)	g	2000 (E2)	1000 (E2)	1000 (E2)	
Application range (temperature)		With "isoCAL" function: +5 +40 °C Without "isoCAL" function: +15 +25 °C			
Display result					
(depending on the set filter level)		0.1 - 0.4			
Weighing pan size (W × D)	mm	140 × 140			
Weighing chamber height (draft shield DE)	mm	172	<u> </u>		
Protection		Protected against du	ust and water		

^{*} DIR = Directive 90/384/EEC on non-automatic weighing instruments used within the European Economic Area

^{** =} Typical min. initial weighing according to USP (Unites States Pharmacopeia), USP31-NF26

Verified r	nodels ง	with EC	Type.	Approval	Certificate:	Precision	balances
------------	----------	---------	-------	----------	--------------	-----------	----------

Models		MSA623S-0CE	MSA623P-0CE	MSA323S-0CE		
Accuracy class*						
For verified models: EG-Type-Approval	Certifica	ate D09-09-015, Type:	MSX			
Scale interval d*	mg	1	1/2/5	1		
Weighing capacity max*	g	620	150/300/620	320		
Calibration value e*	mg	10	10	10		
Min. load min*	mg	20	20	20		
Tare equalization range (subtractive)		≤ 100% from max. weighing capacity				
Application range according to DIR*	g	0.02 - 620	0.02 - 620	0.02 - 320		
Min. initial weight**	g	1.5	1.5	1.5		
Typical stabilization time	S	≤ 0.8	≤ 0.8	≤ 0.8		
Typical measurement time	S	≤ 1	≤ 1	≤ 1		
Application range (temperature)		With "isoCAL" function	on: +5 to +40 °C Without '	'isoCAL" function: +10 to +30 °C		
Display result						
(depending on the set filter level)		0.1 - 0.4				
Weighing pan size (W × D)	mm	140 × 140				
Weighing chamber height (draft shield D	E)) mm	172				
Protection		Protected against de	ust and water	·		

Models		MSA14202S- 0CE	MSA14202P- 0CE	MSA10202S- 0CE	MSA8202S- 0CE
Accuracy class*		I	I	I	
For verified models: EC Type-Approval	Certificat	e D09-09-015, Ty	pe: MSX		
Scale interval d*	g	0.01	0.01/0.02/0.05	0.01	0.01
Weighing capacity max*	g	14,200	3500/7000/ 14,200	10,200	8200
Calibration value e*	g	0.1	0.1	0.1	0.1
Min. load min*	g	1	1	1	0.5
Tare equalization range (subtractive)	re equalization range (subtractive) ≤ 100% from r				
Application range according to DIR*	g	1 - 14,200	1 - 14,200	1 - 10,200	0.5 - 8200
Min. initial weight**	g	15	15	12	12
Typical stabilization time	S	≤ 1	≤ 1	≤ 1	≤ 1
Typical measurement time	S	≤ 1.5	≤ 1.5	≤ 1.5	≤ 1.5
Application range (temperature): With "isoCAL" function Without "isoCAL" function		+5 to +40 °C +15 to +25 °C	+5 to +40 °C +15 to +25 °C	+5 to +40 °C +15 to +25 °C	+5 to +40 °C +10 to +30 °C
Display result (depending on the set filter level)		0.1 - 0.4			
Weighing pan size (W \times D)	mm	206 × 206			
Protection		Protected agains	st dust and water		

^{*} DIR = Directive 90/384/EEC on non-automatic weighing instruments used within the European Economic Area
** = Typical min. initial weighing according to USP (Unites States Pharmacopeia), USP31-NF26

Model-specific Data

Verified models with EC Type Approval Certificate: Precision balances

Models		MSA6202S- OCE	MSA6202P- 0CE	MSA5202S- 0CE	MSA4202S- 0CE
Accuracy class*				I	I
For verified models: EC Type-Approval	Certifica	te D09-09-015, Typ	e: MSX		
Scale interval d*	g	0.01	0.01/0.02/0.05	0.01	0.01
Weighing capacity max*	g	6200	1500/3000/6200	5200	4200
Calibration value e*	g	0.1	0.1	0.1	0.1
Min. load min*	g	0.5	0.5	1	0.5
Tare equalization range (subtractive)		≤ 100% from ma	x. weighing capacity		
Application range according to DIR*	g	0.5 - 6200	0.5 - 6200	1 - 5200	05 - 4200
Min. initial weight**	g	12	12	10	12
Typical stabilization time	S	≤ 1	≤ 1	≤ 0.8	≤ 0.8
Typical measurement time	S	≤ 1.5	≤ 1.5	≤ 1	≤ 1
Application range (temperature): With "isoCAL" function Without "isoCAL" function		+5 to +40 °C +10 to +30 °C	+5 to +40 °C +10 to +30 °C		+5 to +40 °C +10 to +30 °C
Display result (depending on the set filter level)		0.1 - 0.4			
Weighing pan size (W × D)	mm	206 × 206			
Protection		Protected agains	t dust and water		

Models		MSA2202S- 0CE	MSA1202S- 0CE	MSA12201S- 0CE	MSA8201S- OCE	MSA5201S- 0CE
Accuracy class*		II		II	I	II
For verified models: EC Type-Approval	Certificat	e D09-09-015, Ty	pe: MSX			
Scale interval d*	mg	10	10	100	100	100
Weighing capacity max*	g	2200	1200	12,200	8200	5200
Calibration value e*	g	0.1	0.1	1	1	1
Min. load min*	g	0.5	0.5	5	5	5
Tare equalization range (subtractive)		≤ 100% from m	ax. weighing ca	pacity		
Application range according to DIR*	g	0.5 - 2200	0.5 - 1200	5 - 12,200	5 - 8200	5 - 5200
Min. initial weight**	g	12	12	100	100	100
Typical stabilization time	S	≤ 0.8	≤ 0.8	≤ 0.8	≤ 0.8	≤ 0.8
Typical measurement time	S	≤ 1	≤ 1	≤ 1	≤ 1	≤ 1
External standard calibration value (min. accuracy class)	kg	2 (F1)	1 (F1)	10 (F1)	5 (F2)	5 (E2)
Application range (temperature)		With "isoCAL" fu	nction: +5 to +4	O °C Without "	isoCAL" function:	+10 to +30 °C
Display result						
(depending on the set filter level)		0.1 - 0.4				
Weighing pan dimensions (W + D)	mm	206×206				
Protection		Protected agains	st dust and wat	er		


^{*} DIR = Directive 90/384/EEC on non-automatic weighing instruments used within the European Economic Area


** = Typical min. initial weighing according to USP (Unites States Pharmacopeia), USP31-NF26


Balance Dimensions


Microbalances

All dimensions are given in millimeters

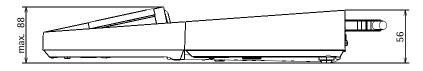

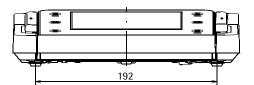


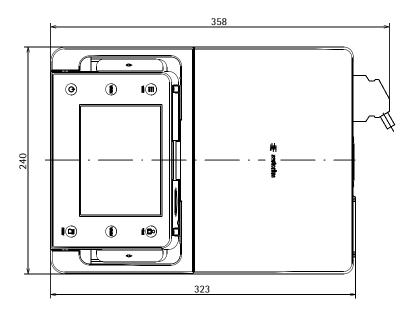

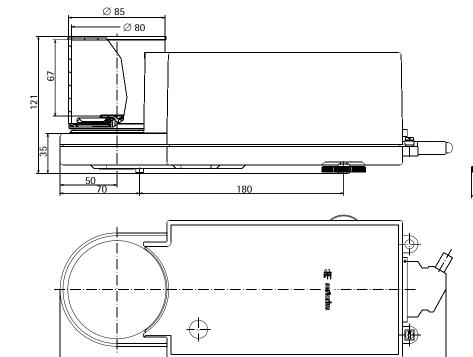


246

340

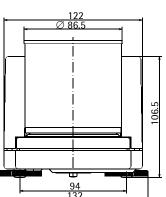

Filter-Microbalances

All dimensions are given in millimeters

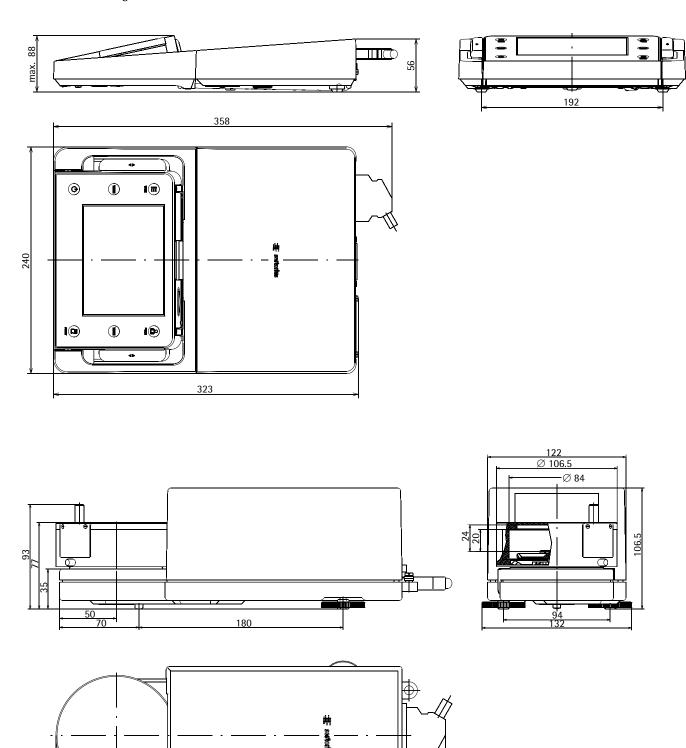


246

344


Semi-microbalances with Motorized Draft Shield


All dimensions are given in millimeters

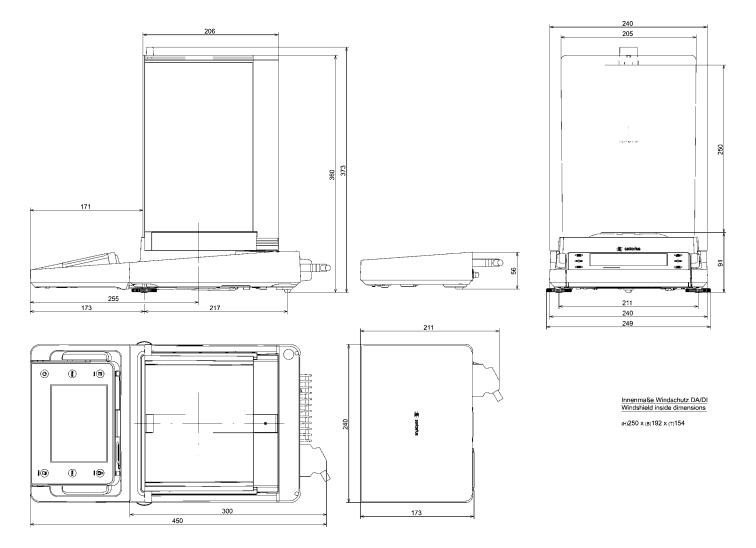


Semi-microbalances and Analytical Balances with Manual DU Draft Shield

All dimensions are given in millimeters

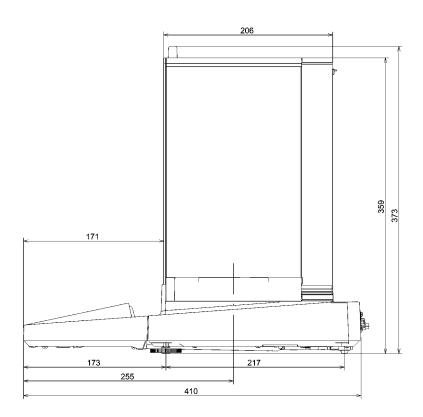

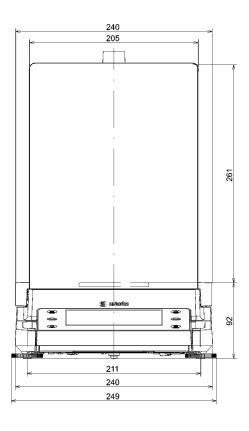


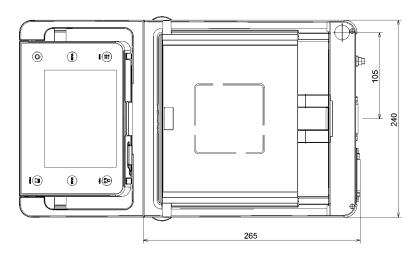



Windshield inside dimensions
(H)261 x (B)193 x (T)191

Precision Balances with a Readability of 1 mg and Manual DE Draft Shield

All dimensions are given in millimeters

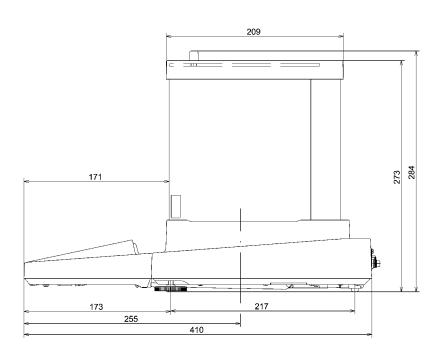

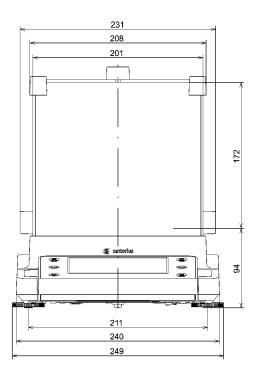


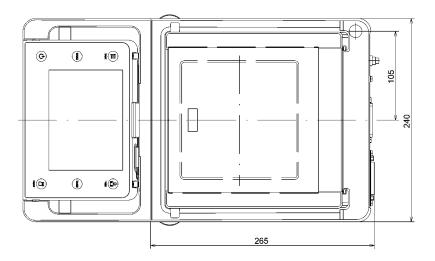



Windshield inside dimensions

(H)172 x (B)193 x (T)191

Precision Balances with a Readability of 1 mg and Framed DR Draft Shield All dimensions are given in millimeters

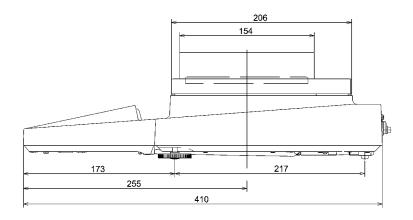


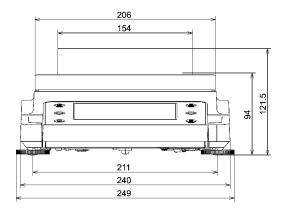


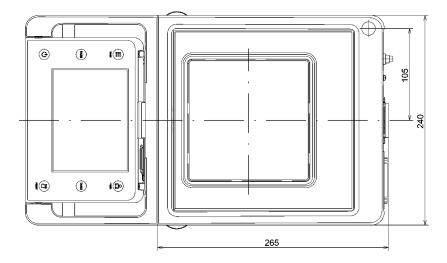

Precision Balances with no Draft Shield

All dimensions are given in millimeters

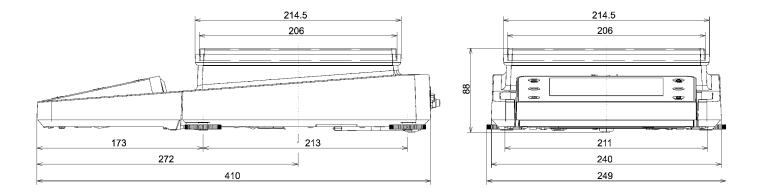

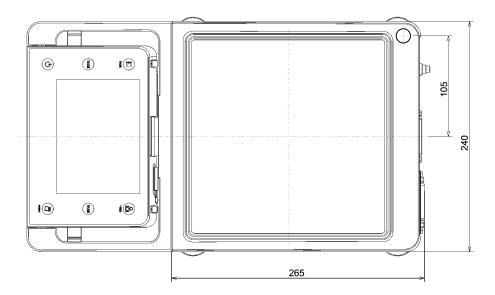

Accessories

Cubis® Optional Accessories

D : 4		^	
Printers	and	Commu	าเดอร์เกท

Printers and Communication	
Verifiable data printer for connection to RS-232, 25-pin accessory interface	YDP10-0CE
Verifiable data printer with <i>Bluetooth</i> * data transmission (with YD001MS-B or IB option only)	YDP10BT-0CI
Color ribbon for YDP10-0CE and YDP10BT-0CE	6906918
Paper rolls for printer YDP10-0CE; 5 rolls 50 m each	6906937
Bluetooth® data interface for wireless connection of data printer YDP10BT	YD001MS-B
RS232C data interface, 9-pin including PS/2 for connecting a PC or keyboard	YD001MS-P
RS232C data interface, 25-pin for connection of Cubis® accessories	YD001MS-R
Display cable 3 m for Cubis® MSA and MSU models, for separate setup of display and weighing unit (Installation by Sartorius Service or in factory [order VF4016])	YCC01-MSD3
Cable 3 m between weighing module and electronics module for Cubis $^{\circ}$ models with 0.01 mg 0.001 mg 0.0001 mg readability	YCC01-MSM3
Installation display cable 3 m for Cubis® models, for separate setup of display and weighing unit	VF4016
RS232C connection cable to connect PC with 9-pin COM interface, length 1.5 m	7357314
SartoCollect software for data communication between balance and PC	YSC02
Sartorius OPC Server for connecting all Sartorius Cubis® balances Requires 32-bit Microsoft Windows 2000 or XP with current service packs. (free download of a 30-day trial version from the Sartorius website) - Initial license - Each additional license within an order	62890PC 62890PC-L
Displays and Input Output Elements	
MSA control unit with color TFT graphic display and touch screen	YAC01MSA
MSE display unit with backlit LC display and tactile keys	YAC01MSE
MSU control unit with backlit b w graphic display and tactile navigation keys	YAC01MSU
Barcode reader with connection cable, 120 mm reading range	YBR03PS2
Foot switch for printing, taring, or using function keys, selection via menu, incl. T connector	YFS01
Infrared sensor for touch-free activation of functions (e.g., draft shield control)	YHS01MS
Hand switch for printing, taring, or using function keys, selection via menu, incl. T connector	YHS02
Foot switch for the draft shield OPEN CLOSED functions (only in combination with DA and DI draft shield), taring and printing	YPE01RC
Additional display, LCD, figure size 13 mm, backlit	YRD03Z
3-segment control display, red – green – red, for plus minus measurements, incl. T connector	YRD11Z

189

Pipette calibration kit (hardware) for models with 0.1 mg and 0.01 mg readability

Consists of moisture trap and all required adapters	
Pipette calibration kit (hardware) for microbalance weighing modules 6.6S and 3.6P	VF988
Consists of moisture trap and all required adapters	
Pipette Tracker pipette calibration software. Software and user manual in English only.	YCP04-PT
Pipette Tracker Pro pipette calibration software, for use in regulated areas, networkable and validatable, according to the 21 CFR Part 11 regulations. Software and user manual in English only.	YCP04-PTPro
Documentation basis for validation (IQ, OQ) of Pipette Tracker PRO version. All documents are in English only.	YCP04-VTK
Filter Weighing and Antistatic Accessories	
Antistatic weighing pan, diameter 130 mm, for weighing modules with a readability of 0.1 mg or 0.01 mg	YWP01MS
Filter weighing pan d 75 mm, for ultramicrobalance or microbalance models (weighing modules 6.6S, 2.7S; only together with DF draft shield)	VF2562
Filter weighing pan d 90 mm, for ultramicrobalance or microbalance models (weighing modules 6.6S, 2.7S; only together with DF draft shield)	VF2880
(weighing modules 6.65, 2.75, only together with Dr draft shield)	
lonization blower to eliminate electrostatic charges on sample containers and samples	YIB01-DR

YCP04MS

Special Applications

Density determination kit for solids and liquids for weighing modules with a readability < 1 mg	YDK01MS
Density determination kit for solids and liquids for weighing modules with a readability of 1 mg	YDK02MS
Q-Grip, flexible holder for weigh-in containers and filters up to 120 mm diameter	YFH01MS
(replaces the original weighing pan, for Cubis® models with 0.01 and 0.1 mg readability)	

Q-Grid grid weighing pan for Cubis® models with a readability of 10 mg or 100 mg for weighing in laboratory hoods, YWP03MS safety weighing cabinets or workbenches (reduced wind attack surface of the weighing pan; replaces the standard weighing pan)

Weighing Tables

Weighing table made from synthetic stone, with vibration dampening	YWT03
Wall console	YWT04
Weighing table made from wood with synthetic stone for precise, reliable measurements	YWT09

Weighing Accessories

Weighing scoop made from chrome nickel steel , 90 + 32 \times 8 mm	641214
Aluminum weighing scoop, 4.5 mg (250 pieces) for ultramicrobalance and microbalance models	6565-250
Aluminum weighing scoop, 52 mg (50 pieces) for ultramicrobalance and microbalance models	6566-50
Support arm for 10/100 mg precision weighing modules for raising the control units MSE, MSU, MSA	YDH01MS

The brand name and logo for *Bluetooth*® wireless technology are are the property of Bluetooth SIG Inc. The use of this brand name and trademark by Sartorius is under license. Other brand names and trademarks are the property of their respective owners.

Declarations of Conformity

Weighing instruments for use in legal metrology: Council Directive 2009/23/EC "Non-automatic Weighing Instruments"

This Directive regulates the determination of mass in legal metrology. For the respective Declaration of Conformity for weighing instruments that have been verified by SARTORIUS for use as legal measuring instruments and that have an EC Type-Approval Certificate, see next page.

This Directive also regulates EC verification by the manufacturer, provided that an EC Type Approval Certificate has been issued and the manufacturer has been accredited by a notified body registered at the Commission of the European Community for performing such verification.

The legal basis for Sartorius to perform the EC verification is EC Directive No. 2009/23/EC for non-automatic weighing instruments. This Council Directive has that has been in effect since January 1, 1993 in the Internal Market. The further legal basis is founded on the Sartorius Quality Management System issued by the Metrology Department of the Regional -Administration Office of Lower Saxony, Germany ("Niedersächsisches Landesverwaltungsamt - Eichwesen") on February 15, 199.

"Installation" Service in Germany

Our "Installation" service package provides a range of important services that guarantee your satisfaction with our work:

- Setup
- Operation
- Inspection
- Instruction

If the initial installation of the weighing instrument is to be carried out by Sartorius, please request this service from a customer service employee.

Re-verification in Germany

The validity of the verification ends when the next but one calendar year has elapsed. When the weighing instrument is used for the control of filling quantities according to the regulation on prepackaging, the verification ends when the next calendar year has elapsed. Re-verification should be requested in good time from the local Weights and Measures office. As appropriate, please observe all statutory amendments.

Subsequent Verifications within the European Countries

The expiration date of the verification depends on the national regulations of the country in which the weighing instrument is used. For information on verification and legal regulations currently applicable in your country, and to obtain the names of the persons to contact, please contact your local Sartorius office, dealer or service center.

Further information concerning verification can be obtained from our customer service centers.





Sartorius Weighing Technology GmbH Weender Landstr. 94 – 108 D-37075 Goettingen, Germany

erklärt, dass das Betriebsmittel declares that the equipment

Geräteart: Device type: Elektronische Halbmikro-, Analysen- und Präzisionswaage

Electronic Semi-micro, Analytical and Precision Balance

Baureihe | Type series:

MSA....- MSE....- MSU....-..

in der von uns in Verkehr gebrachten Ausführung mit den grundlegenden Anforderungen der folgenden Europäischen Richtlinien übereinstimmt:

in the form as delivered complies with the basic requirements of the following European Directives:

Richtlinie 2004/108/EG Directive 2004/108/EC Elektromagnetische Verträglichkeit

Electromagnetic compatibility

Richtlinie 2006/95/EG

Elektrische Betriebsmittel zur Verwendung innerhalb bestimmter

Spannungsgrenzen

Directive 2006/95/EC

Electrical equipment designed for use within certain voltage limits

Richtlinie 1999/5/EG

Funkanlagen und Telekommunikationsendeinrichtungen und die

gegenseitige Anerkennung ihrer Konformität

Directive 1999/5/EC

Radio equipment and telecommunications terminal equipment and the mutual

recognition of their conformity

Das Gerät erfüllt die anwendbaren Anforderungen der in Anhang aufgeführten harmonisierten Europäischen Normen.

The apparatus meets the applicable requirements of the harmonized European Standards listed in Annex.

Jahr der Anbringung der CE-Kennzeichnung / Year of attachment of CE marking: 11

Sartorius Weighing Technology GmbH Goettingen, 2011-06-30

Dr. Reinhard Baumfalk Vice President R&D

Dr. Dieter Klausgrete

Leitung International Certification Management Head of International Certification Management

Diese Erklärung bescheinigt die Übereinstimmung mit den genannten EG-Richtlinien, ist jedoch keine Zusicherung von Eigenschaften. Bei einer mit uns nicht abgestimmten Änderung des Produktes verliert diese Erklärung ihre Gültigkeit. Die Sicherheitshinweise der zugehörigen Produktdokumentation sind zu beachten. This declaration certifies conformity with the above mentioned EC Directives, but does not guarantee product attributes. Unauthorised product modifications make this declaration invalid. The safety information in the associated product documentation must be observed.

SWT11CE001

34785-000-58

SOP-3.RD-045-fo2



EG-Konformitätserklärung EC Declaration of Conformity

Anhang / Annex

Liste der angewendeten harmonisierten Europäischen Normen List of the applied harmonized European Standards

Richtlinie 2004/108/EG / Directive 2004/108/EC

EN 61326-1:2006 Elektrische Mess-, Steuer-, Regel- und Laborgeräte - EMV- Anforderungen - Teil 1: Allgemeine Anforderungen (IEC 61326-1:2005)

Electrical equipment for measurement, control and laboratory use - EMC requirements - Part 1: General requirements (IEC 61326-1:2005)

Richtlinie 2006/95/EG / Directive 2006/95/EC

EN 61010-1:2001 Sicherheitsbestimmungen für elektrische Mess-, Steuer-, Regel- und Laborgeräte - Teil 1: Allgemeine Anforderungen (IEC 61010-1:2001)

Safety requirements for electrical equipment for measurement, control, and laboratory use - Part 1: General requirements (IEC 61010-1:2001)

3. Richtlinie 1999/5/EG / Directive 1999/5/EC

Nur für Geräte mit Bluetooth Datenausgang Typ YBT03

Only for devices with Bluetooth data output interface type YBT03

EN 301 489-1 V1.8.1 Elektromagnetische Verträglichkeit und Funkspektrumangelegenheiten (ERM) — Elektromagnetische Verträglichkeit (EMV) für Funkeinrichtungen und -dienste — Teil 1: Gemeinsame technische Anforderungen Electromagnetic compatibility and Radio spectrum Matters (ERM) — ElectroMagnetic Compatibility (EMC) standard for radio equipment and services — Part 1: Common technical requirements

EN 301 489-17 V1.3.2 Elektromagnetische Verträglichkeit und Funkspektrumangelegenheiten (ERM) — Elektromagnetische Verträglichkeit für Funkeinrichtungen und dienste — Teil 17: Spezifische Bedingungen für Breitbandübertragungssysteme im 2,4 GHz Band, Einrichtungen in lokalen Hochleistungs-Funk-

netzen (RLAN) im 5 GHz Band und Breitband-Datenübertragungssysteme

im 5,8 GHz Band

Electromagnetic compatibility and Radio spectrum Matters (ERM) — ElectroMagnetic Compatibility (EMC) standard for radio equipment — Part 17: Specific conditions for 2,4 GHz wideband transmission systems, 5 GHz high performance RLAN equipment and 5,8 GHz Broadband Data Transmitting Systems

EN 300 328 V1.7.1 Elektromagnetische Verträglichkeit und Funkspektrumangelegenheiten (ERM) — Breitband-Übertragungssysteme — Datenübertragungsgeräte, die im 2,4-GHz-ISM-Band arbeiten und Breitband-Modulationstechniken verwenden — Harmonisierte EN, die wesentliche Anforderungen nach

Artikel 3 Absatz 2 der R&TTE-Richtlinie enthält

Electromagnetic compatibility and Radio spectrum Matters (ERM) — Wideband transmission systems — Data transmission equipment operating in the 2,4 GHz ISM band and using wide band modulation techniques — Harmonized EN covering essential requirements under Article 3(2) of the REITE Directive

.

SWT11CE001 34785-000-58 SOP-3.RD-045-fo2





Sartorius Weighing Technology GmbH Weender Landstr. 94 – 108 37075 Goettingen, Germany

erklärt, dass das Betriebsmittel declares that the equipment

Geräteart: Device type: Elektronische Halbmikro-, Analysen- und Präzisionswaage

Electronic Semi-micro, Analytical and Precision Balance

Baureihe / Type series:

MSA....- MSE....- MSU....- MSU....-

übereinstimmt mit den Regelungen der Europäischen Richtlinie (in der heute gültigen Fassung): complies with the regulations of the European Directive (in the today valid version):

Richtlinie 2002/95/EG

Zur Beschränkung der Verwendung bestimmter gefährlicher Stoffe

in Elektro- und Elektronikgeräten

Directive 2002/95/EC

on the restriction of the use of certain hazardous substances in electrical

and electronic equipment

sofern das Betriebsmittel gekennzeichnet ist mit: provided that the equipment is marked with:



Sartorius Weighing Technology GmbH Goettingen, 2011-06-24

Dr. Reinhard Baumfalk

Leitung, Entwicklung, Mechatronik Vice President, R&D, Mechatronics Dr. Dieter Klausgrete

Leitung International Certification Management, Mechatronik Head of International Certification Management, Mechatronics

Diese Erklärung bescheinigt die Übereinstimmung mit der genannten EG-Richtlinie, ist jedoch keine Zusicherung von Eigenschaften. Bei einer mit uns nicht abgestimmten Änderung des Produktes verliert diese Erklärung ihre Gültigkeit

This declaration certifies conformity with the above mentioned EC Directive, but does not guarantee product attributes. Unauthorised product modifications make this declaration invalid.

SWT11RoHS001

34785-000-58

SOP-3.RD-045-fo3

C E Declaration of Type Conformity to Directive No. 2009/23/EC

This declaration is valid for non-automatic electromechanical weighing instruments for use in legal metrology. These weighing instruments accepted for legal metrological verification have an EC Type-Approval Certificate. The model(s) concerned is (are) listed below along with the respective type, accuracy class, and EC Type-Approval Certificate number:

Model	Weighing instrument type	Accuracy class	EC type-approval certificate no.
MSCE	MSX	0	D09-09-015
GBB,,CE	MSX	0	D09-09-015
MSCE	MSX		D09-09-015

Type weighing module: SA EA, SB EA, SC EA, SD EE, SE EA, SF EA, SB EB, SC EB, SC EC, SG EE, SH EE Type indicating and operator terminal: YACO1MSA, YACO1MSE, YACO1MSU

SARTORIUS Weighing Technology GmbH * declares that its weighing instrument types comply with the requirements of the Council Directive on non-automatic weighing instruments, no. 2009/23/EC of 23 April 2009; the associated European Standard "Metrological aspects of nonautomatic weighing instruments," No. EN 45501; the most recently amended versions of the national laws and decrees concerning legal metrology and verification in the Member States of the European Union, the EU, and the Signatories of the Agreement on the European Economic Area, which have adopted this Council Directive into their national laws; and with the requirements stipulated on the Type-Approval Certificate for verification. This Declaration of Type Conformity is valid only if the ID label on the weighing instrument has the CE mark of conformity and the green metrology sticker with the letter

"M" stamped on it (the two-digit number in large print stands for the year in which the mark was affixed):

Example (date/year and number of the notified body may vary):

C € 11 m

If these marks are not on the ID label, this Declaration of Type Conformity is not valid. Validity can be obtained, for example, by submitting the weighing instrument for final processing by an authorized representative of SARTORIUS Weighing Technology GmbH *. The period of validity of this Declaration of Type Conformity shall expire upon any tampering with, repair or modification of this weighing instrument or, in some Member States, on the date of expiration. This declaration applies only to the weighing instrument without peripheral devices. The operator of this weighing instrument shall be responsible for obtaining an authorized renewal of the verification, such as subsequent or periodic verification, of the weighing instrument for use as a legal measuring instrument.

Sartorius Weighing Technology GmbH * 37070 Goettingen, Germany Signed in Goettingen on 05 August 2011

Dr. Reinhard Baumfalk (Vice President R&D)

J. Rehwald (Head of the Production Department

Mechatronics / Weighing Technology Division)

*) Sartorius Weighing Technology GmbH is the legal successor of Sartorius AG

LOP-3,225_an2e_2005.06.09.doc P106ep03.doc

Physikalisch-Technische Bundesanstalt



Braunschweig und Berlin



EG-Bauartzulassung

EC type-approval certificate

Zulassungsinhaber:

Sartorius Weighing Technology GmbH

Issued to:

Weender Landstr. 94-108

37075 Göttingen

Rechtsbezug: In accordance with: Richtlinie 2009/23/EG vom 23. April 2009 über nichtselbsttätige

Waagen (ABI. L 122 S. 6), umgesetzt durch die Eichordnung vom 12. August 1988, zuletzt geändert am 06.06.2011 (BGBI. I S. 1035). Directive 2009/23/EC of 23 April 2009 on non-automatic weighing instruments (OJ L 122 p. 6), implemented by the Verification Ordinance of 12 August 1988, last

amended on 06.06.2011 (Federal Law Gazette I, p. 1035).

Bauart:

Nichtselbsttätige elektromechanische Präzisionswaage Non-

automatic electromechanical high accuracy weighing instrument

In respect of: Typ/Type:

Max 14200 g, e = 1...200 mg, n ≤ 520000

(II) Max 12200 g, e = 0,01...2 g, n ≤ 82000

Option: Mehrteilungswaage / multi-interval instrument Mehrbereichwaage / multi-range instrument

Zulassungsnummer:

D09-09-015 3. Revision

Approval number:

18.06.2019

Gültig bis: Valid until

Anzahl der Seiten:

Number of pages.

PTB-1.12-4052066

Geschäftszeichen: Reference No.:

Benannte Stelle:

0102

Notified Body:

Ort, Ausstellungsdatum:

Braunschweig, 07.07.2011

Date of issue:

Zertifizierer:

Certifier: Im Auftrag

By order

Timo Schwabe





Bewerter: Evaluator: Im Auftrag By order


Dipl.-Ing. M. Link

Die Hauptmerkmale, Zulassungsbedingungen und Auflagen sind in der Anlage enthalten, die Bestandteil der Revision der EG-Bauartzulassung ist. Hinweise und eine Rechtsbehelfsbelehrung befinden sich auf der ersten Seite der Anlage

The principal characteristics, approval conditions and special conditions, if any, are set out in the Annex which forms an integral part. of this Revision of the EC type-approval certificate. For notes and information on legal remedies, see first page of the Annex.

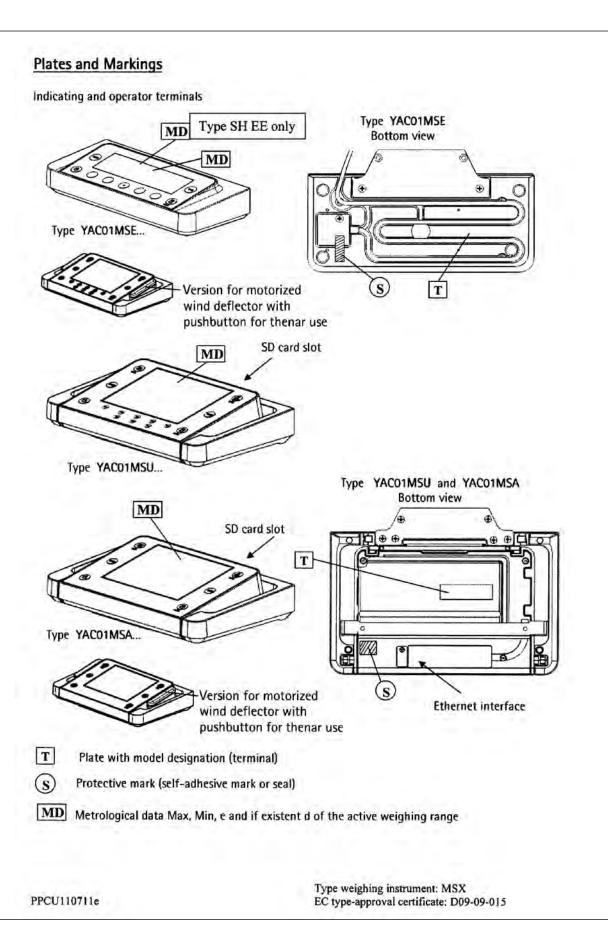

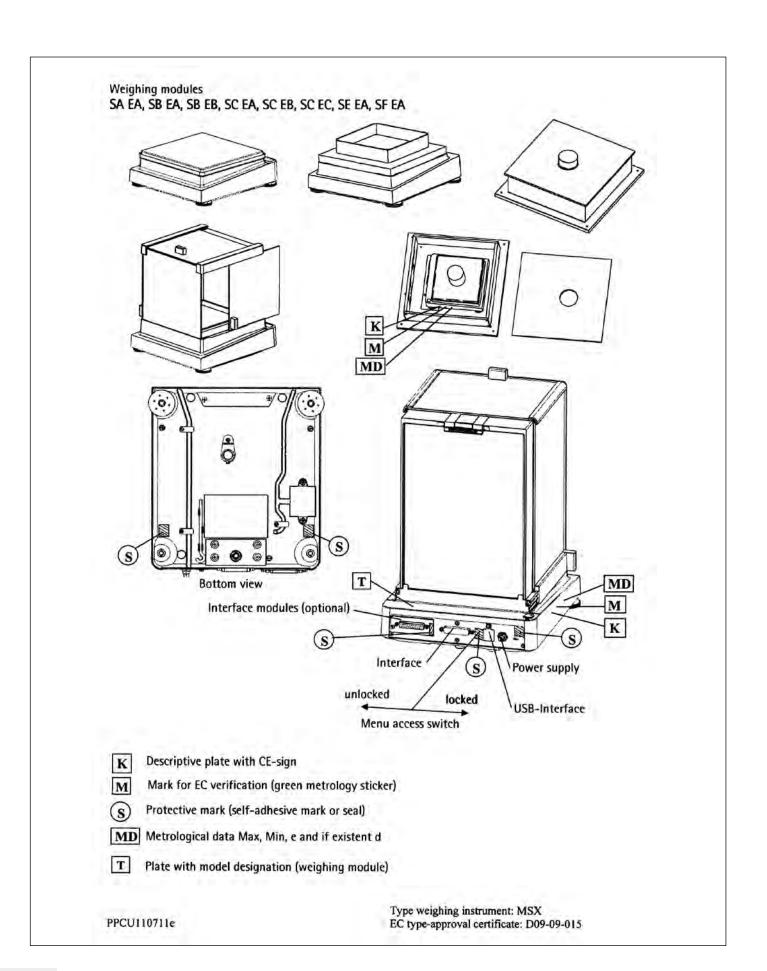

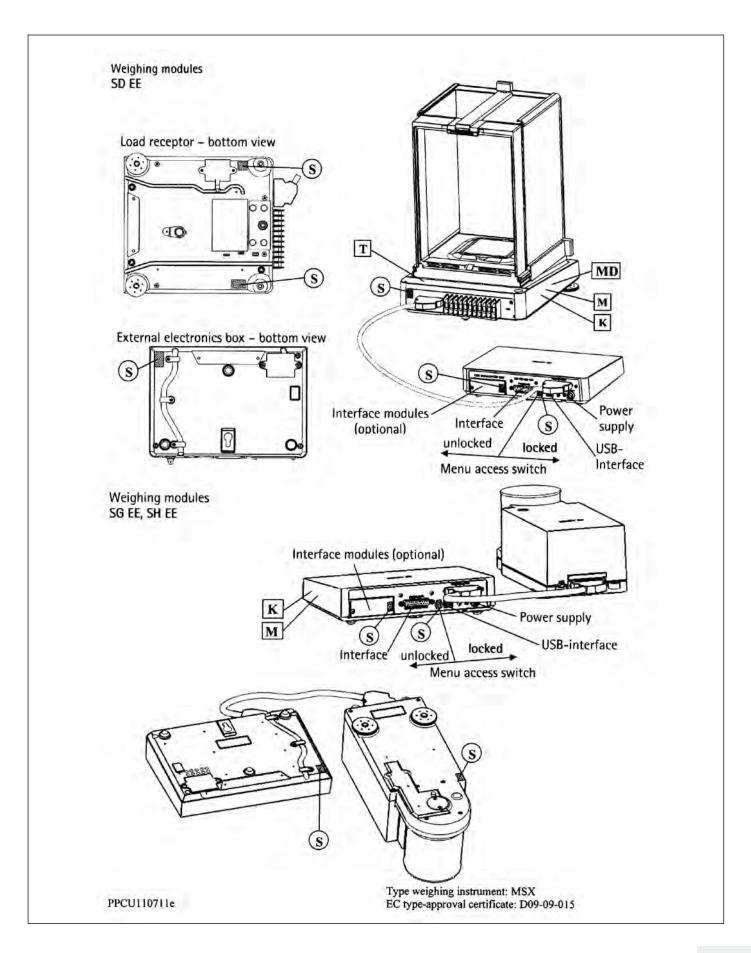

Type MSX

Indicating and operator terminals: YAC01MSE, YAC01MSA, YAC01MSU Weighing modules: SA EA, SB EA, SB EB, SC EA, SC EB, SC EC, SD EE, SG EE, SH EE

Example of descriptive plate on a weighing instrument already verified **K**



Sartorius Weighing Technology GmbH,	D09-09-015 MSX (SC EC) Max	12200 g	
Goettingen, Germany	+10°C/+30°C	Min	5 g	
C€09 M	+ 5°C./+40°C iso-CAL	e =	1 g	
CC03 0111 1VI	II 1111444	d =	0,1 g	

Example of plate with model designation (weighing module) T



Example of plate with model designation (terminal) T

Sartorius Weighing Technology GmbH, Goettingen, Germany

Nicht unter Last verbinden oder trennen! YAC01MSE

Do not connect or disconnect when energized!

PPCU110711e

Type weighing instrument: MSX EC type-approval certificate: D09-09-015

Sartorius Weighing Technology GmbH Weender Landstrasse 94–108 37075 Goettingen, Germany

Phone +49.551.308.0 Fax +49.551.308.3289 www.sartorius.com

Copyright by Sartorius, Goettingen, Germany. All rights reserved. No part of this publication may be reprinted or translated in any form or by any means without the prior written permission of Sartorius.

The status of the information, specifications and illustrations in this manual is indicated by the date given below.

Sartorius reserves the right to make changes to the technology, features, specifications and design of the equipment without notice.

Status:

November 2011, Sartorius Weighing Technology GmbH, Goettingen, Germany