


Sample Banking Management System for iSBS

# SBMS 2.0

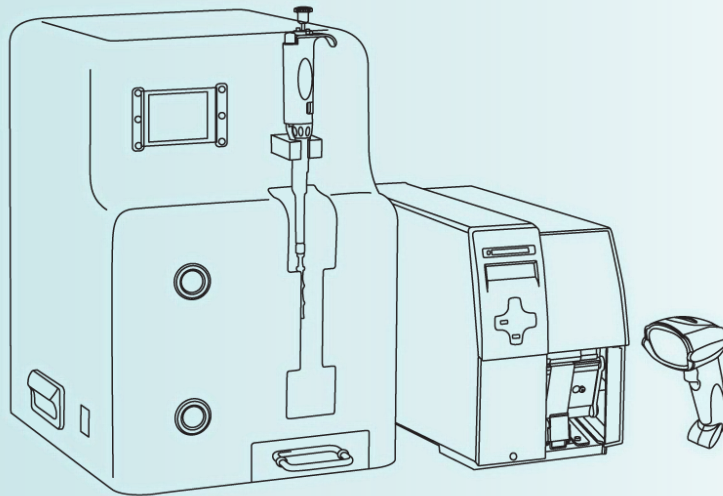

## Installation & User Manual

Version 2.1

**MicroDigital Co., Ltd.**  
#A-101, Korea Bio Park B/D, 694-1, SamPyung-Dong,  
BunDang-Gu, SungNam-Si, GyungGi-Do, 463-400, Korea  
Tel : +82-31-628-2200 / Fax : +82-31-628-2222  
[info@md-best.com](mailto:info@md-best.com) / [www.MD-Best.com](http://www.MD-Best.com)





## Table of Contents

Chapters	Sections	Pages
1. Hardware Preparation & IP Address / Gateway Setup	1) Hardware Package	4
	2) IP Address and Gateway Setup for Windows XP	5
	3) IP Address and Gateway Setup for Windows 7	7
2. SBMS Software: Installation & ODBC Setup	1) Installation	9
	2) ODBC Setup for Windows XP	12
	3) ODBC Setup for Windows 7	15
3. SBMS Software Setting & Run	1) Run Program	18
	2) Login	18
	3) Master Window (Main Menu)	19
	4) Registration	21
	5) Search	24
	6) Pulling	25
	7) Export to Excel	27
	8) Configuration	28
4. Note		30


### Attention!

Contents of SBMS can be modified without prior notification for performance upgrades.


# 1. Hardware Preparation and IP Address / Gateway Setup

## 1) Hardware Package

### 1.1 iSBS Package

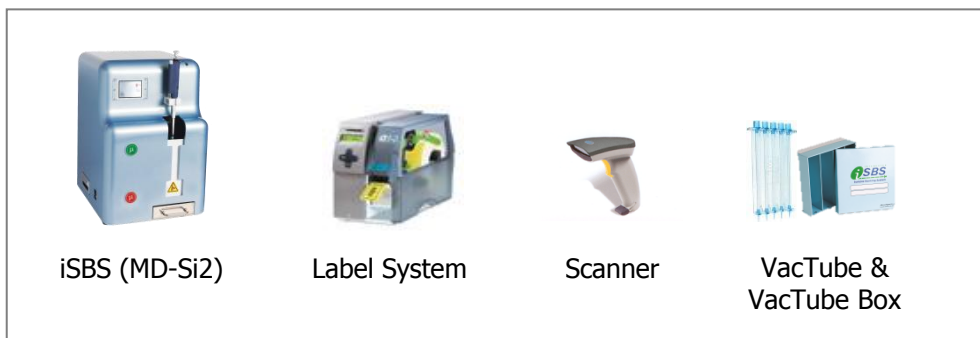

### 1.2 Hardware Setup

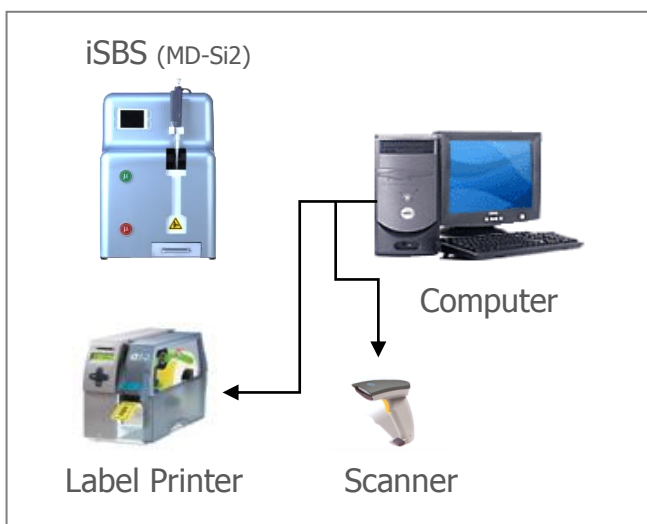


#### **CAUTION**


The computer and the label printer must be connected with Cross LAN Cable, which is included in iSBS Package.

- A. Connect Barcode Printer to your computer via LAN cable (Cross Cable)
- B. Connect Barcode Scanner to Computer via USB port
- C. Power On Barcode Printer
- D. Power On Computer
- E. Set IP Address and Gateway


## 2) IP Address and Gateway Setup for Windows XP

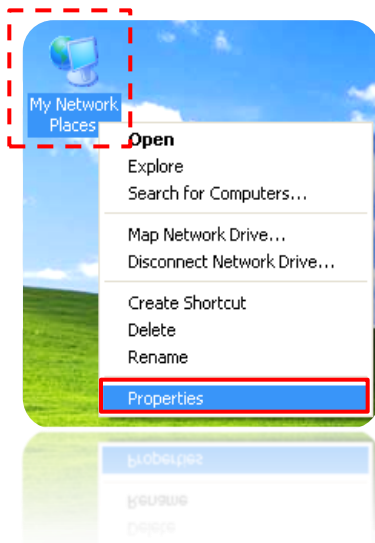


- ① Right click on "My Network Places" from your desktop
- ② Click "Properties" to open the "Network Connections" window

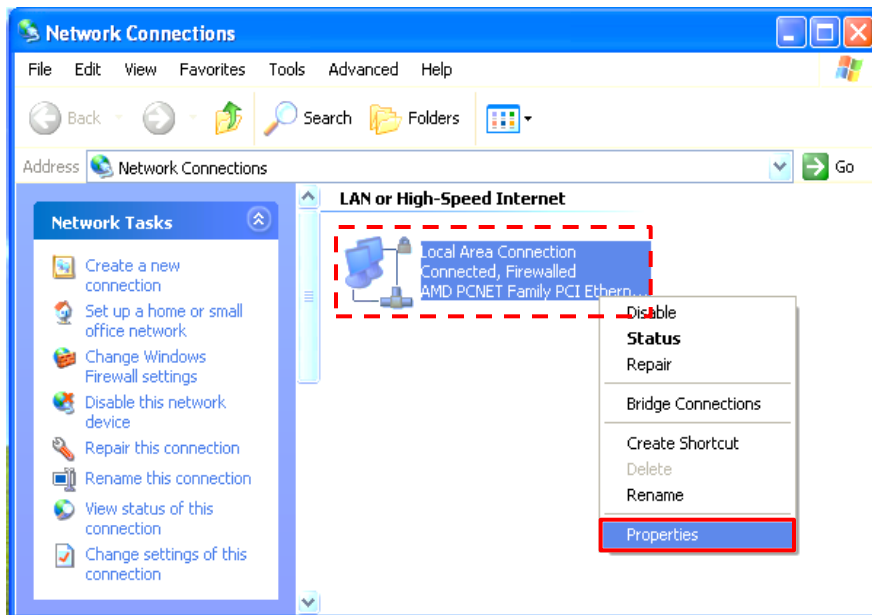

- ③ Right click "Local Area Connection"
- ④ Click "Properties" to open the "Local Area Connection Properties" window

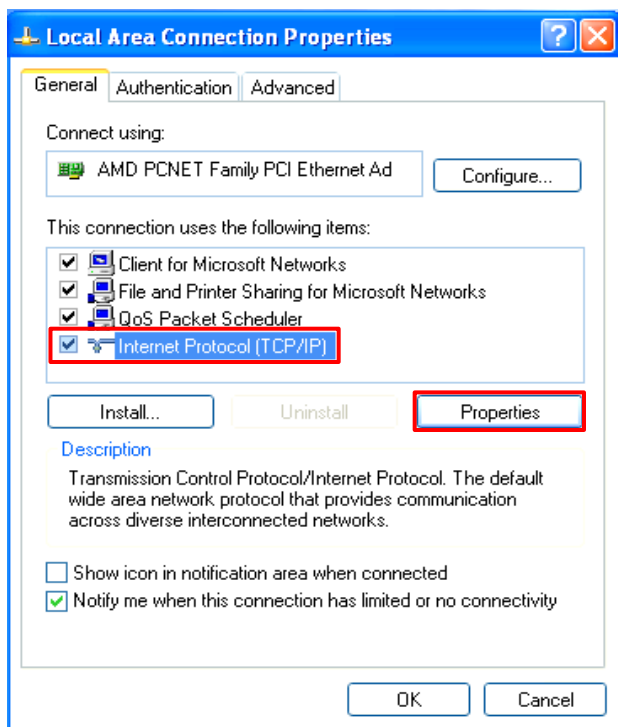

- ⑤ Click "Internet Protocol (TCP/IP)"
- ⑥ Click "Properties" to open "Internet Protocol (TCP/IP) Properties" Window

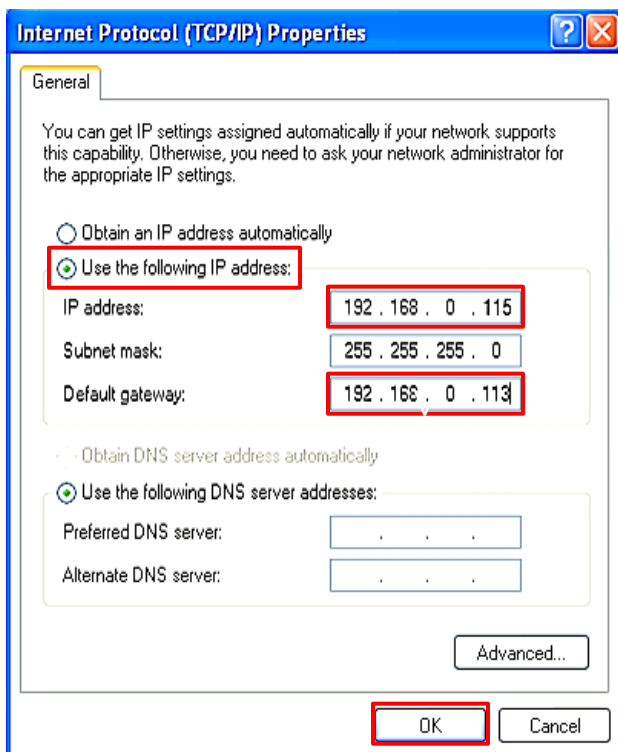


- ⑦ Click and select "Use the following IP address:"
- ⑧ Set "IP address:" to 192.168.0.115
- ⑨ "Subnet mask" values will be automatically appear (Do NOT change!)
- ⑩ Set "Default gateway" to 192.168.0.113 and click "OK" to save and close


### Attention!

The IP address MUST NOT be same as the IP address set in the Label Printer; however Default gateway MUST be same as the Default gateway in the Label Printer.

### 3) IP Address and Gateway Setup for Windows 7

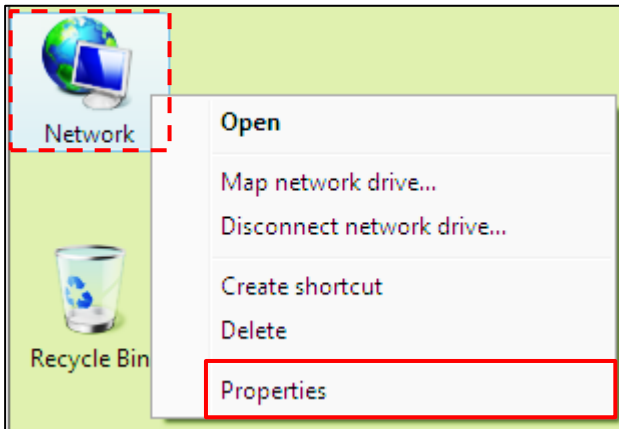


- ① Right click on "Network" from your desktop
- ② Click "Properties" to open up the "Network" window

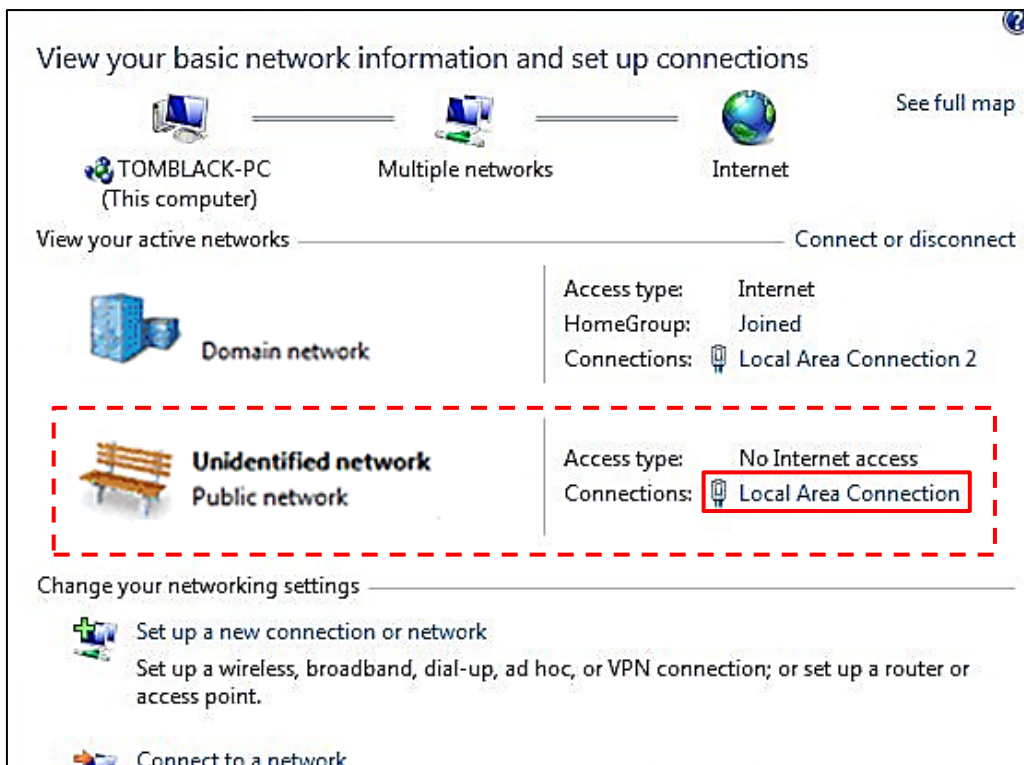


- ③ Click "Local Area Connection" on the right side of "Unknown network" to open Local Area Connection Status window

(Caution: Make sure the label printer is connected to the computer via LAN Cable)

**Attention!** If "Unidentified network" cannot be found, please check the hardware connection between the Label Printer and Computer.

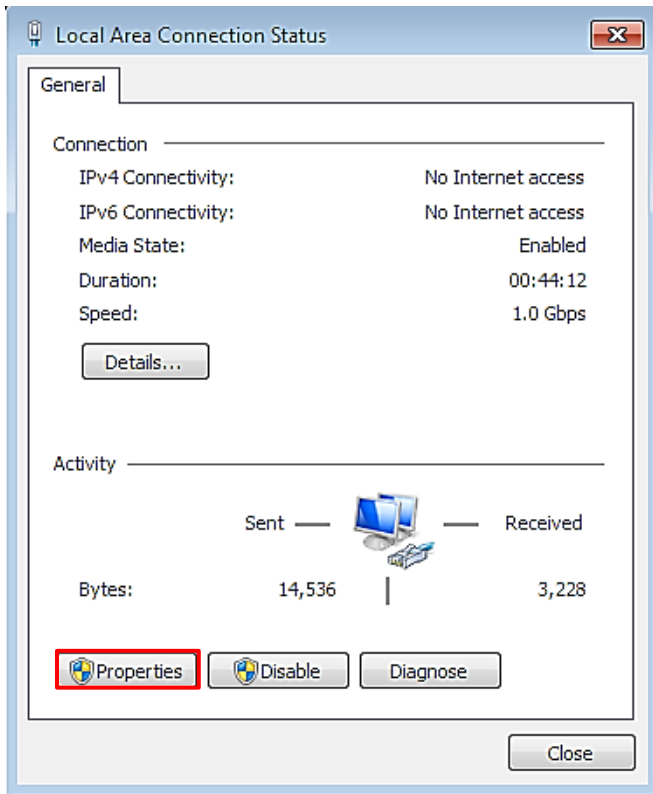


- ④ Click "Properties" in the Local Area Connection Status window

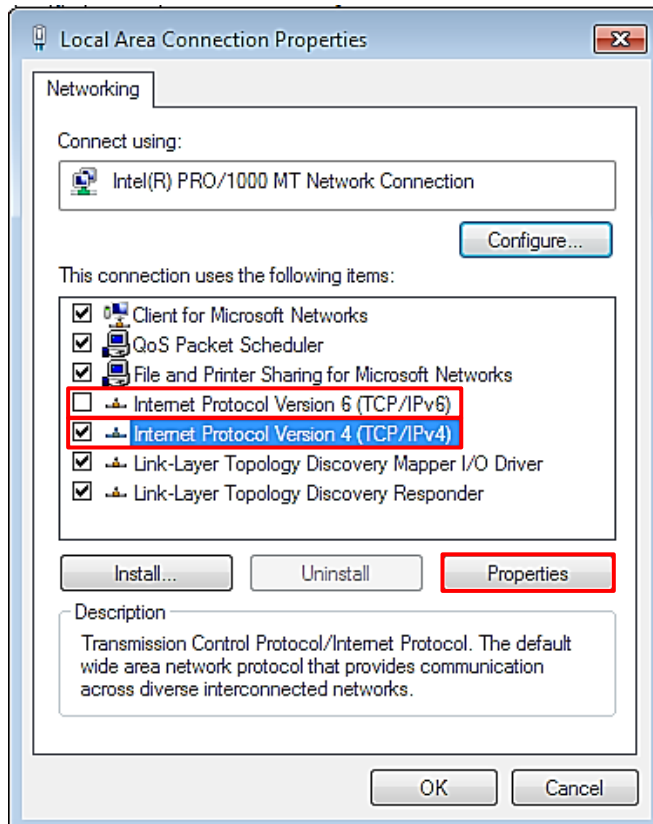

- ⑤ Click and uncheck "Internet Protocol Version 6 (TCP/IPv6)" on the list and make sure "Internet Protocol Version 4 (TCP/IPv4)" is checked
- ⑥ Select "Internet Protocol Version 4 (TCP/IPv4)"
- ⑦ Click "Properties" to open "Internet Protocol Version 4 (TCP/IPv4) Properties" window



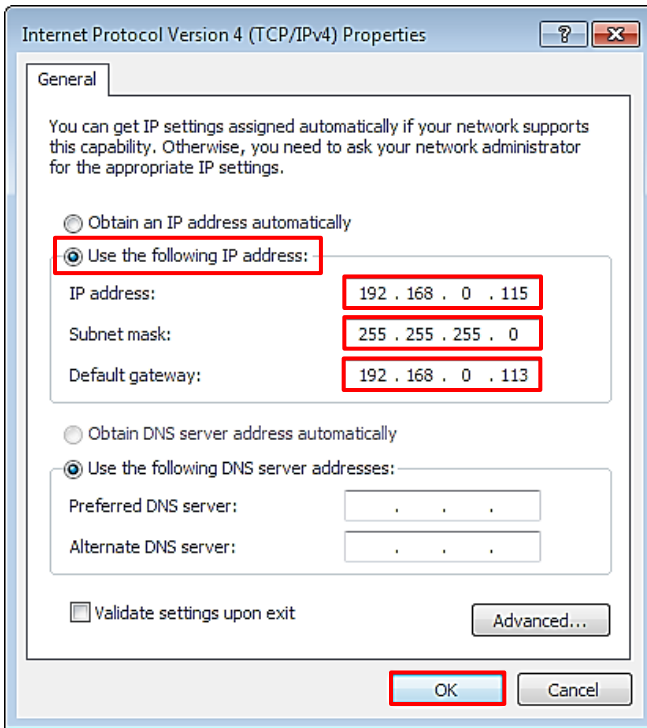

- ⑧ Select "Use the following IP Address : " and set "IP Address" to "192.168.0.115"  
"Subnet Mask" values will automatically appear (Do NOT touch)
- ⑨ Set "Default Gateway" to "192.168.0.113"
- ⑩ Click "OK" and save and close

## 2. SBMS Software: Installation & ODBC Setup

### 1) Installation


#### For Windows XP

- ① Double click on "SBMS (ver2.0)" icon in SBMS CD to start the installation of SBMS Software



#### For Windows 7

- ① Right click on "SBMS (ver2.0)" in the CD
- ② Click on "Run as administrator"

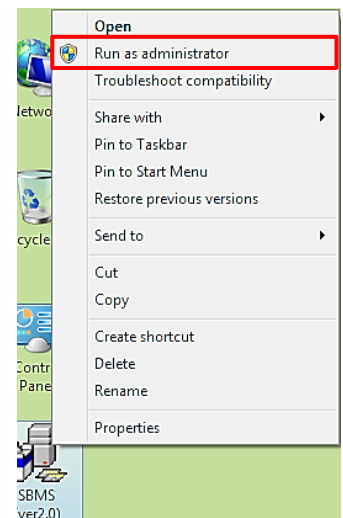




③ Click "Next" to proceed

**Attention!**

It is recommended to exit all other programs before the installation.

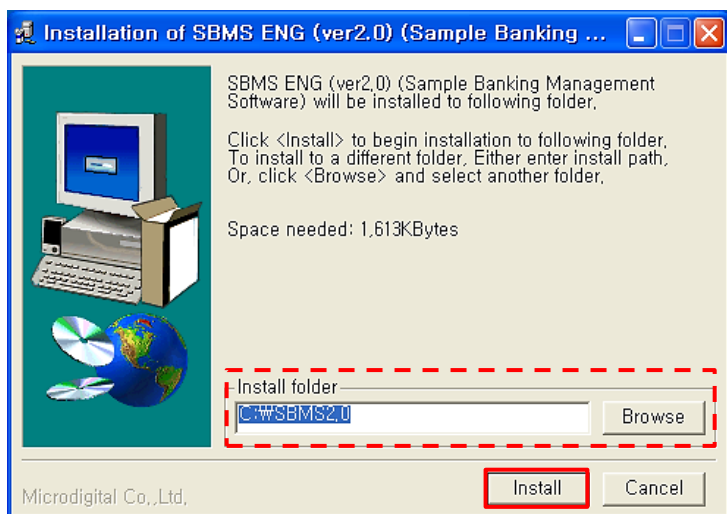

④ Set directories for "Install folder"

⑤ Click "Install" to proceed the installation

**Attention!**

Default Install folder is "C:\ SBMS2.0"






- ④ Click "Ok" to close the window

### Installation complete!

### Attention!

For Windows 7 and VISTA users,  
Please continue to additional steps below

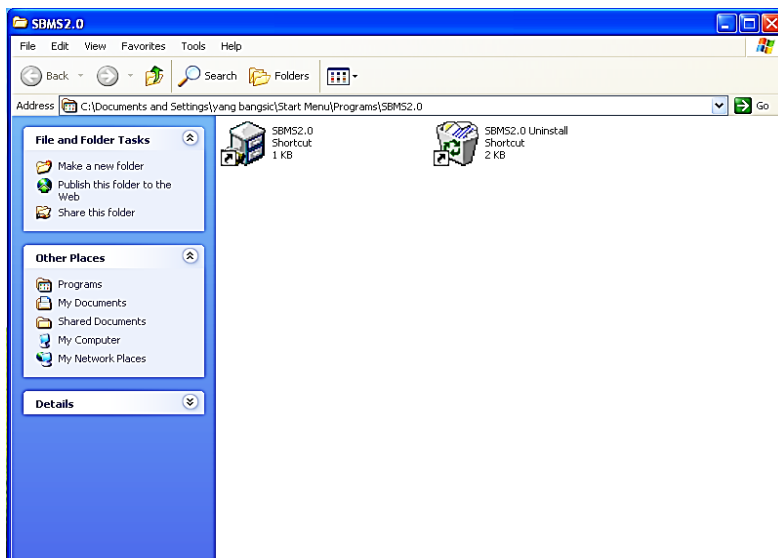


### Attention!

Default Shortcut Location on Start Menu:  
Start → Program → SBMS2.0

### Attention! : Additional steps for Windows 7 and VISTA

- ⑤ Locate LPr.zip file in the provided CD
- ⑥ Extract the LPr.zip file to "SBMS2.0" folder

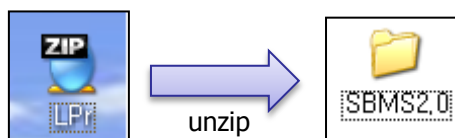


### CAUTION!

Make sure the files compressed in LPr.zip file have transferred directly to "SBMS2.0" folder,  
NOT to any subdirectories in "SBMS2.0" folder.

## 2) ODBC Setup for Windows XP


**Attention!** ODBC needs to be set before execute the program.

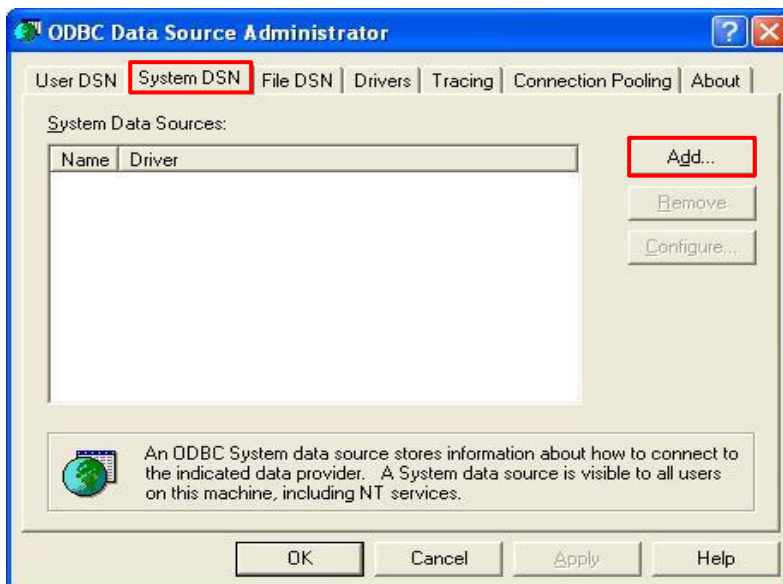

- ① Click "Start → Control → Management Tools → ODBC" to open "ODBC Data Source Administrator" window
- ② Click "System DSN" tab to open
- ③ Click "Add" to open "Create New Data Source" window



- ④ Select "Microsoft Access Driver (\*.mdb)" and click "Finish" to open "ODBC Microsoft Access Setup" window

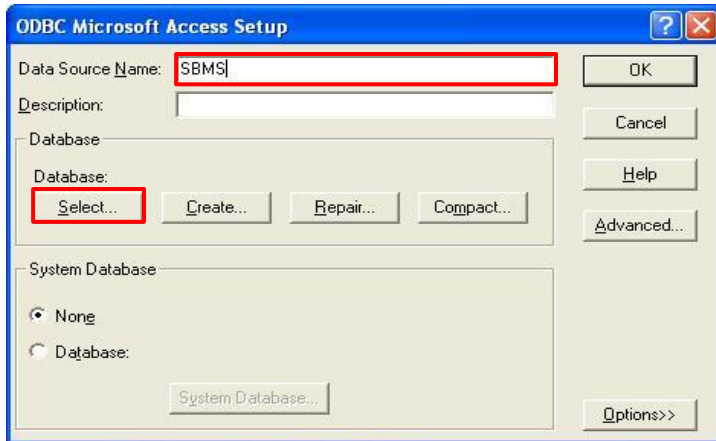


- ⑤ Type "SBMS" in "Data Source Name:"
- ⑥ Click "Select..." to open "Select Database" window




- ⑦ Select "DB\_SBMS.mdb" file in "DATABASE" folder in "SBMS2.0" folder
- ⑧ Click "OK" to return to "ODBC Microsoft Access Setup" window

**Attention!** Default Location for SBMS2.0: C:/SBMS2.0/DATABASE

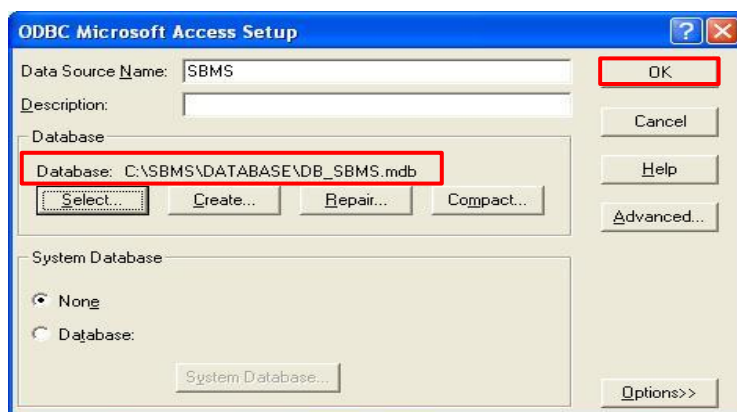

- ⑨ Confirm that DB\_SBMS.mdb file is selected, and click "Ok" to return to "ODBC Date Source Administrator" window

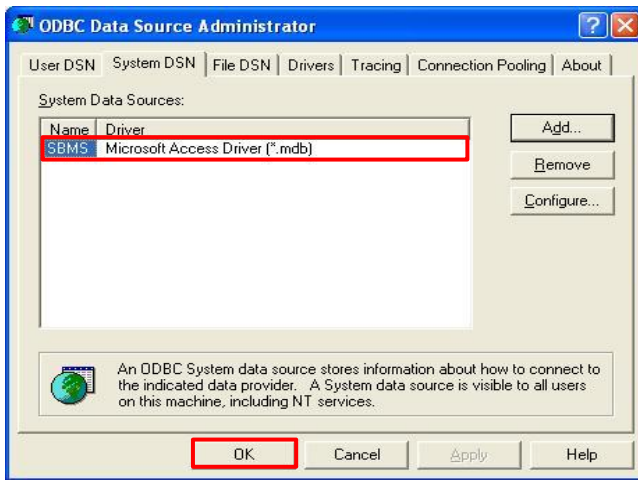


- ⑩ Confirm "SBMS" on the list and click "Ok" to complete the ODBC setup

### 3) ODBC Setup for Windows 7


#### For Windows 7: 32bit

- ① Click "Start" button
- ② Type "ODBC" in search, and press "Enter" to open "ODBC Data Source Administrator" window

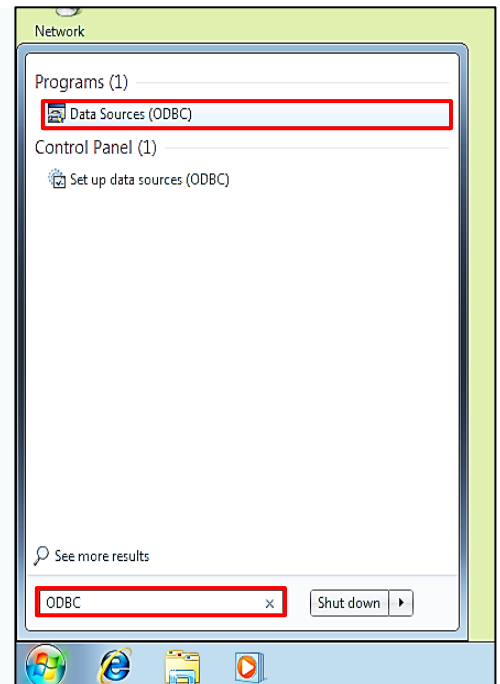

#### For Windows 7: 64bit

- ① Open a "Windows Explorer"
- ② Go to "SysWOW64" folder in "Windows" folder
- ③ Double click "odbcad32" application to open "ODBC Data Source Administrator" window

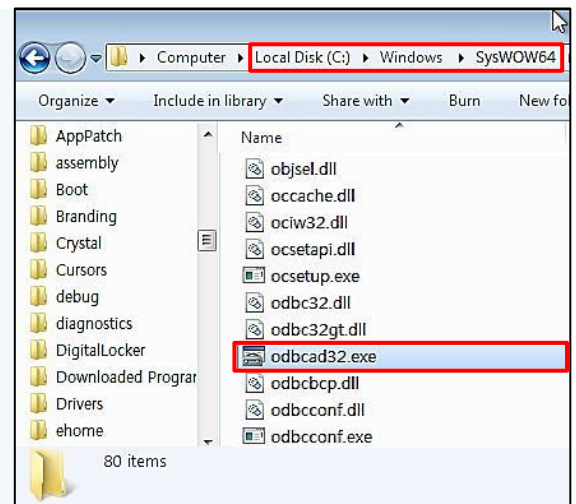


#### CAUTION!


For ODBC setup in Windows: 64bit, users may open "ODBC Data Source Administrator" by following the steps for Windows 7: 32bit above, but you CANNOT add proper contents; therefore make sure to follow proper protocol.

**Attention!**


After the "ODBC Data Source Administrator" window is open, please follow steps below:

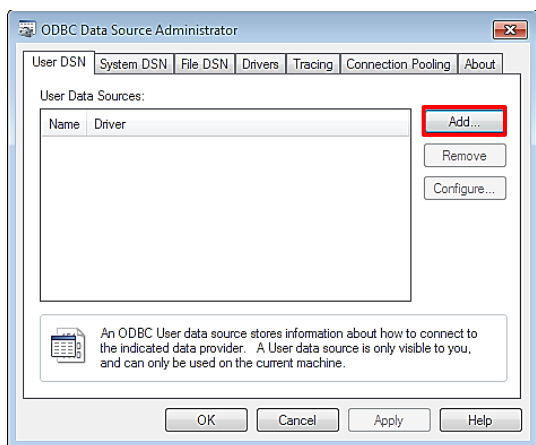

- ① Click "Add" button to open "Create New Data Source" window

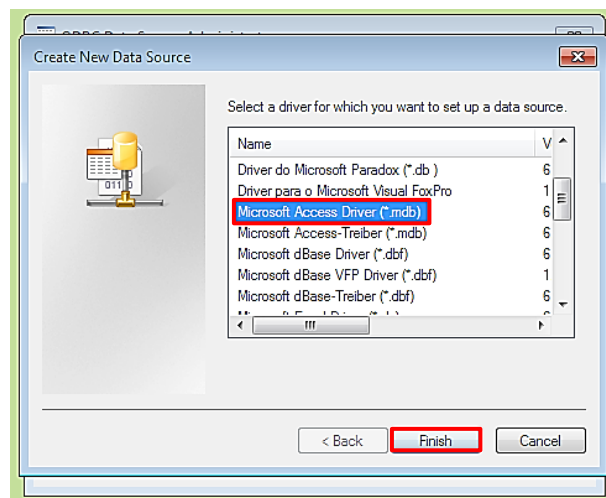


- ② Select "Microsoft Access Driver (\*.mdb)" from the list
- ③ Click "Finish" to proceed to "ODBC Microsoft Access Setup" window

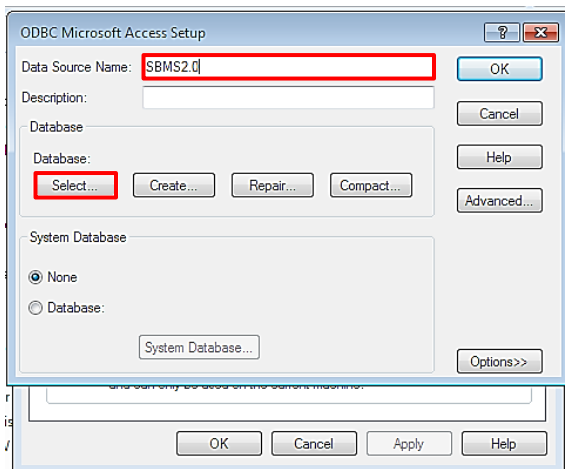

- ④ Type "SBMS2.0" for Data Source Name
- ⑤ Click "Select" to open "Select Database" Windows



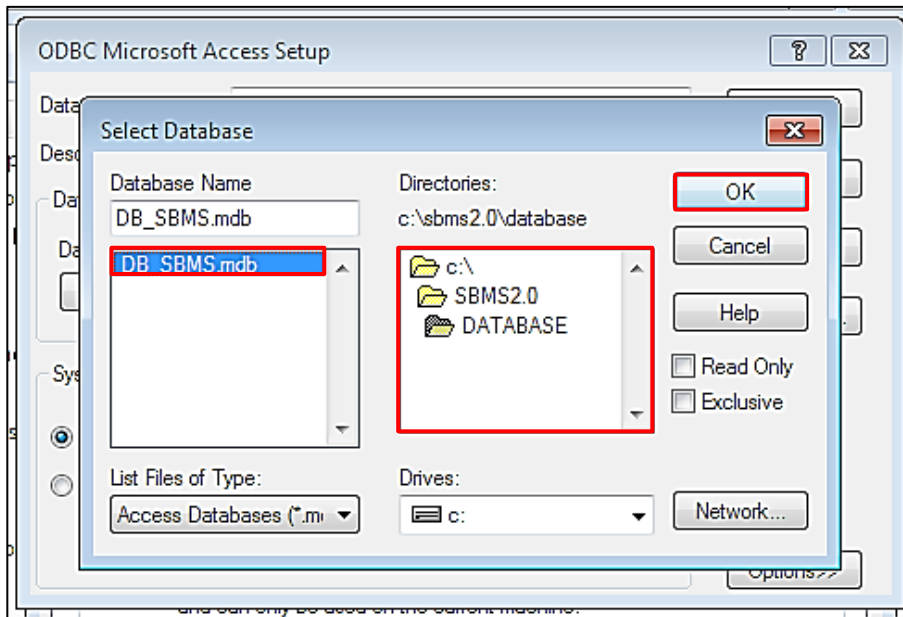

- ⑥ Select "DB\_SBMS.mdb" file in "DATABASE" Folder in "SBMS2.0" Folder  
(Default Location : C:/SBMS2.0/DATABASE)
- ⑦ Close all the windows by clicking "OK"

## 3. SBMS Software Setting & Run

### 1) Run Program

- ① Double clicking “SBMS2.0” icon will run the SBMS software.



#### Attention!

For Windows 7 or VISTA users,


- I. Right click on “SBMS2.0” icon.
- II. Click “Properties” to open “Properties” window for SBMS 2.0 application.
- III. Click “Compatibility” to go to “Compatibility” tab
- IV. Click and check “Run this program as an administrator” under “Privilege Level” Section.

### 2) Login



- First time “Login” requires “User ID” and “Product No.”
- If you do not have them, contact the distributor or manufacturer.
- Password is NOT required for the first time.
- You will set the User ID and Password in the “Config” of SBMS program.


### 3) Master Window (Main Frame)

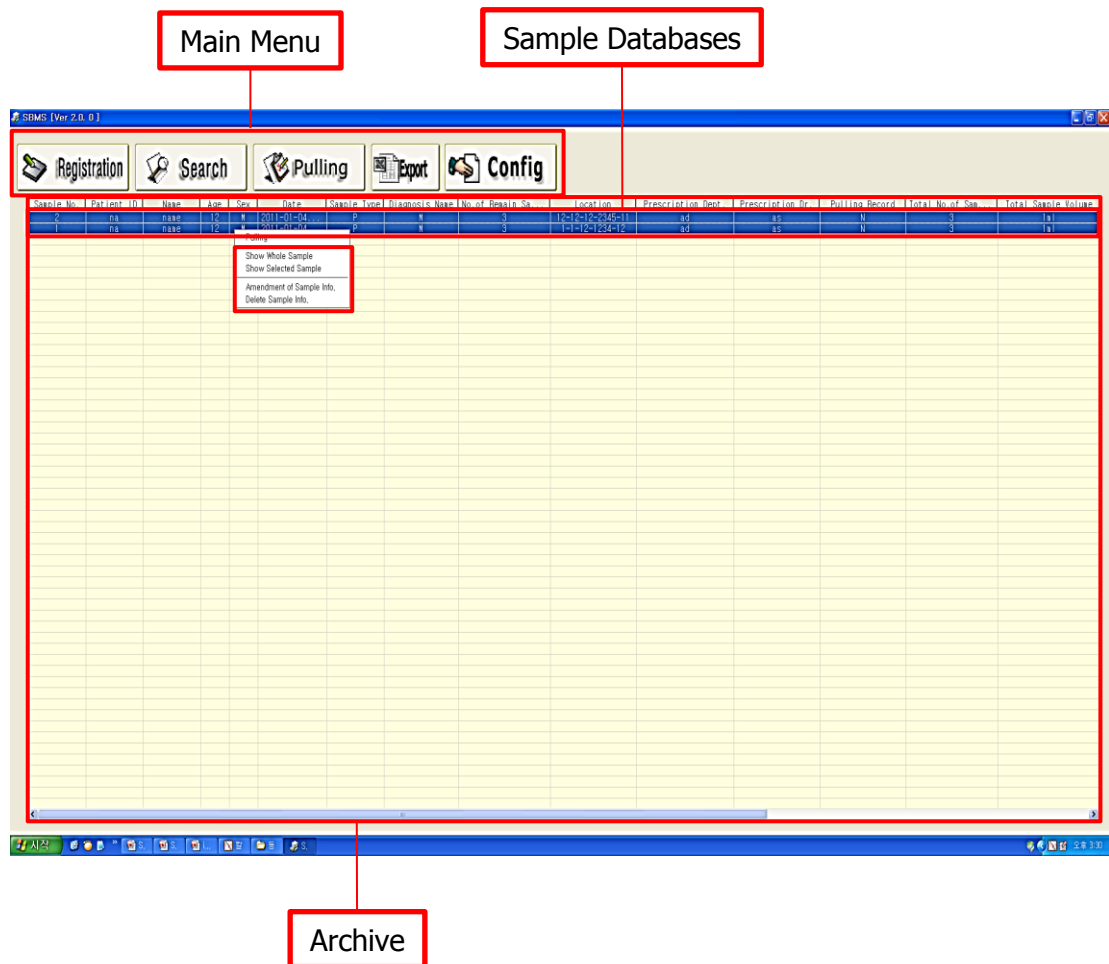

- **Registration** is to register Sample Databases to the Archive.
  - **Search** is to find Sample Databases based on user's search condition.
  - **Pulling** is to update the record of the samples when users pull the samples.
  - **Export** is to export the list of Sample Databases in the Archive to Microsoft Excel.
  - **Config** is to administrate the Archive.
- Only 100 most recently registered Sample Databases are shown in the Archive.
  - Users may select multiple number of Sample Databases by pressing "Ctrl" and select the Databases in the Archive.

## ① Show Whole Samples / Show Selected Samples

- Users can organize the list in the Archive by showing only selected Sample Databases in the Archive by selecting the Sample Databases, right-clicking on the selected Sample Databases, and clicking on “Show Selected Samples”
- To return to the whole list of the Sample Databases in the Archive, right-click on the Archive and select “Show Whole Samples”

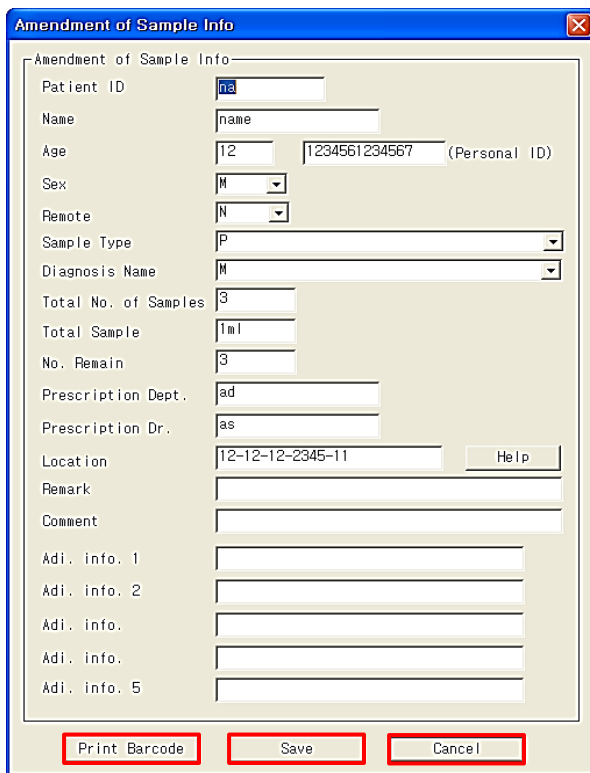


## ② Amendment of Sample Databases

- Users can modify or correct the information in Sample Databases by right-clicking the Sample Database in the Archive and click “Amendment of Sample Info.”
- After the amendment, click “Save” to save the current information in the Database, and click “Print Barcode” to print the barcode for the modified sample information.

- Users may cancel the modification by clicking “Cancel” before save.

**CAUTION!** Users CANNOT print out the barcode without executing save first.

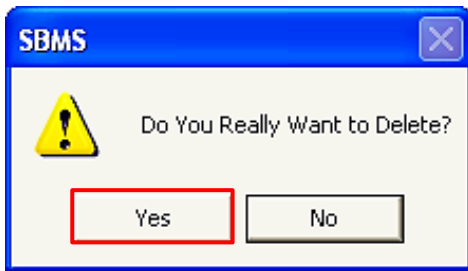


### ③ Deletion of Sample Databases

- Users can delete Sample Databases by right-clicking the Database in the Archive and click “Delete Sample Info.” to open a warning window.  
Click “Yes” to delete the Sample Databases permanently.

## CAUTION!

When the sample databases are deleted, users CANNOT restore the databases.

## 4) Registration

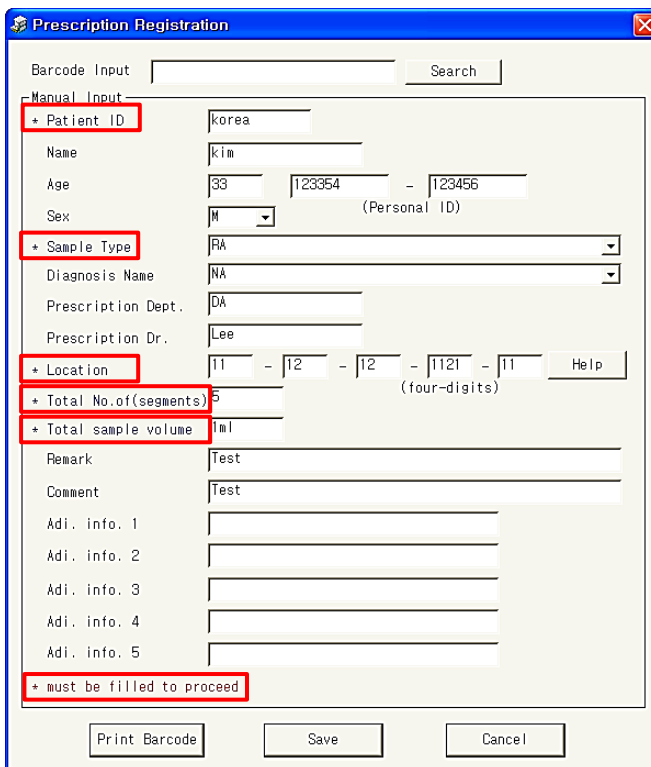

Barcode Input  Search

Manual Input

\* Patient ID

Name

Age   -

Sex  (Personal ID)

\* Sample Type

Diagnosis Name

Prescription Dept.

Prescription Dr.

\* Location  -  -  -  -  Help

\* Total No. of segments  (four-digits)

\* Total sample volume

Remark

Comment

Adi. info. 1

Adi. info. 2

Adi. info. 3


Adi. info. 4

Adi. info. 5

\* must be filled to proceed

### ① Fill in each categories with information of the sample

- Users can register Sample Databases by clicking on “Registration” from Main Menu to open “Prescription Registration” window
- Users can put information of the sample in “Prescription Registration” window.
- Categories with “\*” is required to fill in.
- Users can also search Sample Databases by scanning the barcode from the sample to fill barcode values in “Barcode Input” and clicking on search.

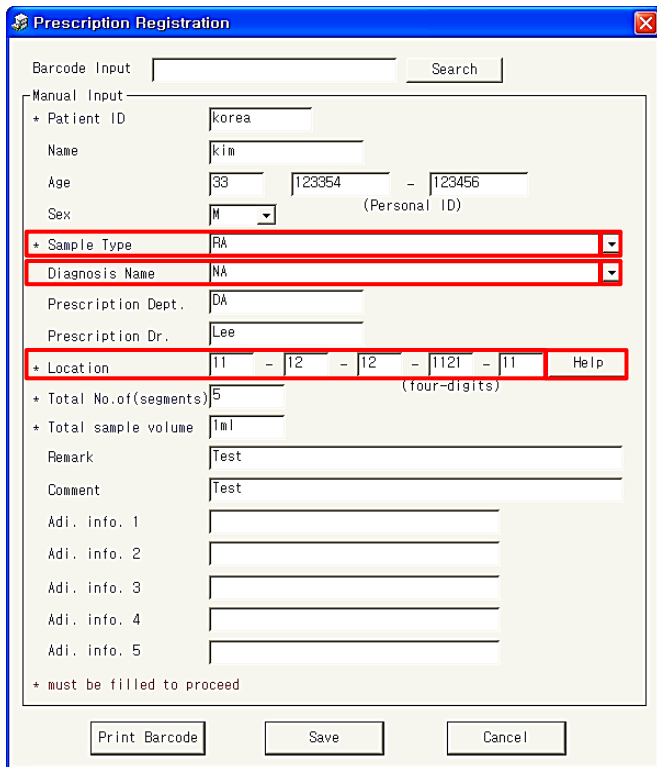


- Users can fill in "Sample Type" manually, however users can also assign Sample Types in Config to select them by scrolling down.
- Users can also assign Diagnosis Names in Config to select them by scrolling down.
- When users click on any boxes for "Location", values will automatically appear, which are the values from the last registration, therefore users can follow the last Sample Database registration.
- If users have selected assigned sample type in "Sample Type" category, values from the 1<sup>st</sup> box in "Location" will be automatically set according to the assigned Sample Type number.


## Attention!

Preregistration (assigning) of Sample Types in Config is recommended.

It is much simpler to manage the "Location" values from the vast amount of Sample Databases in this way.

- Click "Help" to read the description for each box for "Location".
- Values from 1<sup>st</sup> box defines the type of samples.
- Values from 2<sup>nd</sup> box defines the type of deep-freezer. (Deep-freezer ID)
- Values from 3<sup>rd</sup> box defines the location of VacTube Box inside the deep-freezer.
- Values from 4<sup>th</sup> box defines VacTube Box ID.
- Values from 5<sup>th</sup> box defines the location of the sample inside the VacTube Box.

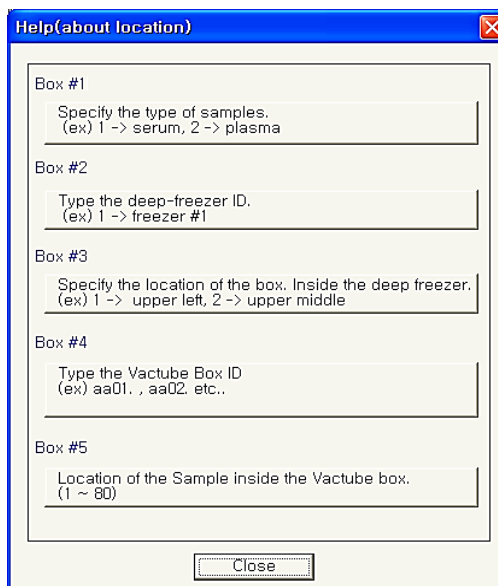


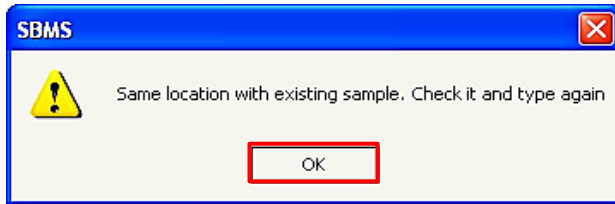

- If the user sets "Location" values identical to other sample which has already been registered, an error occurs.
- In this case, click "OK" to continue the registration process.

## Attention!

Errors occur, if users do not fill in mandatory categories with "\*" Please, make sure to fill in mandatory categories with "\*" before save. Mandatory categories are "Patient ID", "Sample Type", "Location", "Total No.of(segments)", and "Total sample volume"

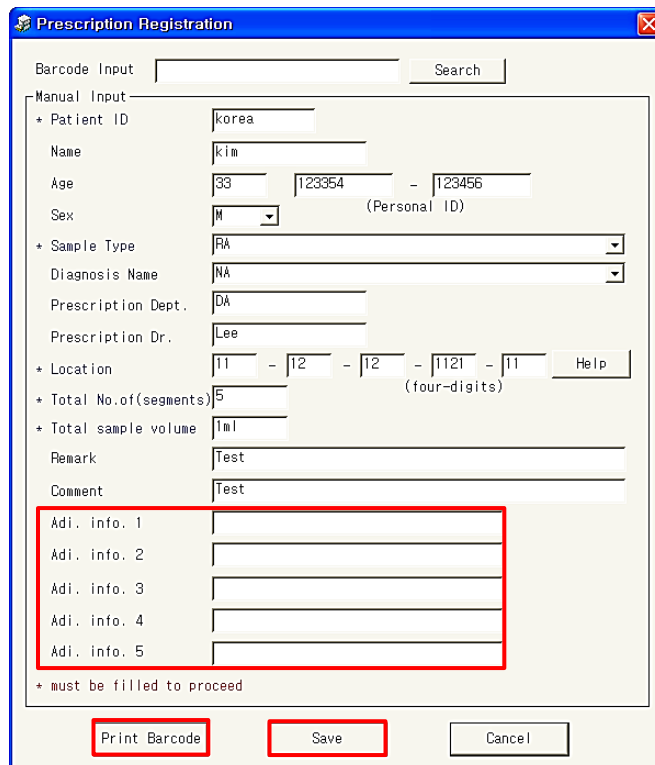

Barcode Input  Search

Manual Input

+ Patient ID

Name

Age   -  (Personal ID)

Sex

+ Sample Type

Diagnosis Name

Prescription Dept.

Prescription Dr.

+ Location  -  -  -  -  Help (four-digits)

+ Total No.of(segments)

+ Total sample volume

Remark

Comment

Adi. info. 1

Adi. info. 2

Adi. info. 3

Adi. info. 4


Adi. info. 5

\* must be filled to proceed

Print Barcode Save Cancel

- For "Adi. Info. #", users can add their own categories by registering categories in Config.
- ② Click "Save" to save current information
  - ③ Click "Print Barcode" to print 2D barcode from Label Printer
  - ④ Attach the printed label to
  - ⑤ Confirm the registered sample database displayed on the uppermost row in the Archive

## 5) Search

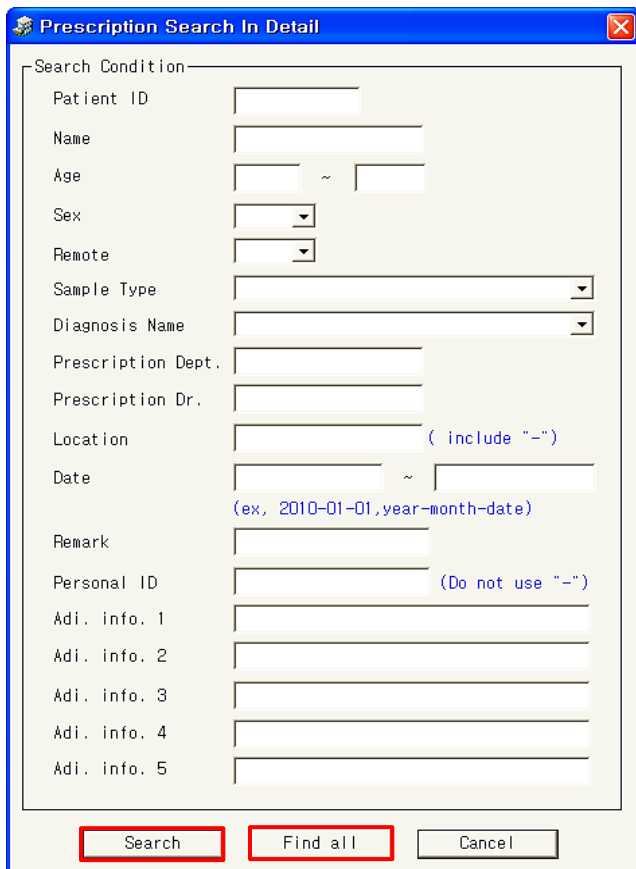

① Click "Search" from Main Menu to open "Prescription Search In Detail" window

② Fill in the categories which will be the condition to search

③ Click "Search" to find matched sample databases based on the categories that are filled in

④ Confirm the matched sample databases

**In order to return to whole list of samples...**

⑤ Click "Search" from Main Menu to open "Prescription Search In Detail" window


⑥ Empty all the categories blank

⑦ Click "Find all" to return to the whole list in the Archive

- In order to search Sample Databases in a certain date, only first blank next to "Date" needs to be filled with the specific date.
- In order to search Sample Databases in a certain age, only first blank next to "Age" needs to be filled with the specific age.
- For all the categories except "Age", "Sex", and "Remote", search system allows users to find any Sample Databases containing the word(s) that the users filled in, within the category. For example, if the user typed "This is" in "Adi. info. 1" category, the user will find any Sample Databases such as Databases containing "This is Sample", and "This is a test" in their "Adi. info. 1", if there are such Sample Databases.



## 6) Pulling

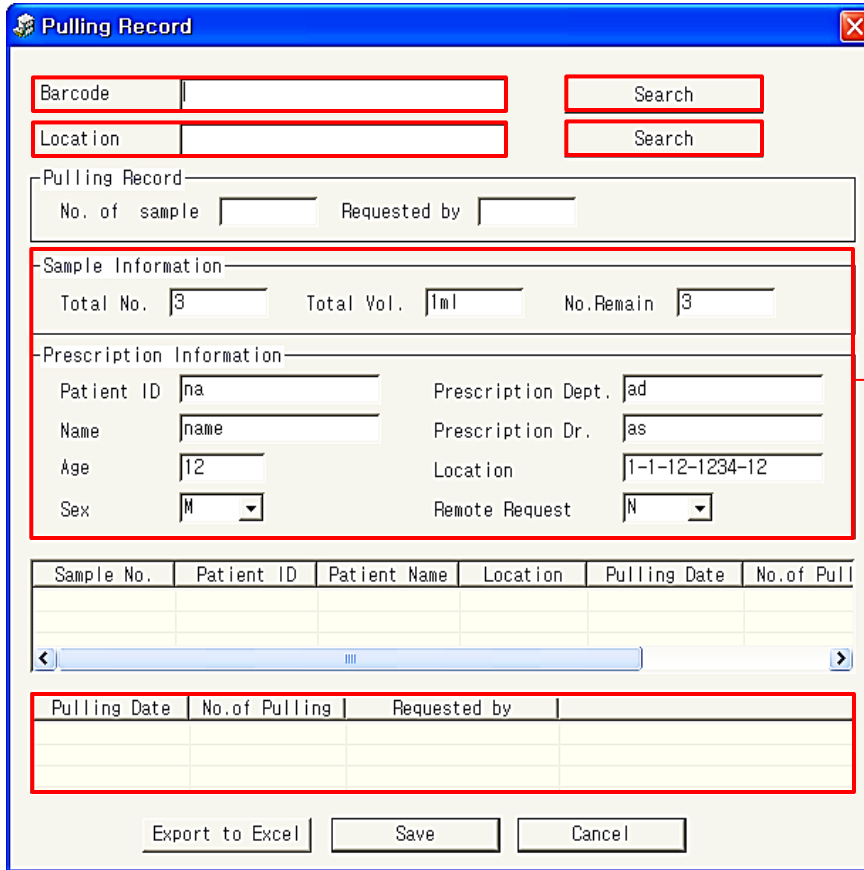

**Pulling Record**

Barcode

Location

-Pulling Record-

No. of sample  Requested by

-Sample Information-

Total No.  Total Vol.  No. Remain

-Prescription Information-

Patient ID  Prescription Dept.

Name  Prescription Dr.

Age  Location

Sex  Remote Request

Sample No.	Patient ID	Patient Name	Location	Pulling Date	No. of Pull

Pulling Date	No. of Pulling	Requested by

Information of the sample

- ① Click "Pulling" from Main Menu to open "Pulling Record" window
- ② Scan the 2-D barcode attached on the VacTube
- ③ Click "Search" to bring out the information of the sample

- In case the user is not able to scan the barcode on the VacTube, user can either manually type in the location code written on the barcode label and click "Search", or select the Database among Sample Databases in the Archive, then right-click, and select "Pulling" to open "Pulling Record" manually.

④ Fill in the categories under "Pulling Record"

⑤ Click "Save"

⑥ Confirm the pulling information

Records of pulling



- Temporary Records of Pulling is available until the user closes the window.
- An error occurs, if there are not enough segments to pull.

- Users can always check the pulling records of samples in "Pulling Record" window.
- In order to amend or delete pulling records in "Pulling Record" window, right-click on the records at the bottom, and select "Delete Sample Info." or "Amendment of Sample Info."
- When the user select "Delete Sample Info.", a confirmation window will be shown.
- Click "Export to Excel" to export the Temporary Records Of Pulling to Microsoft Excel Spreadsheet.



- Select "Yes" to delete the selected pulling records permanently.

H14									
	A	B	C	D	E	F	G	H	I
1	Sample_No	Patient_I	Patient_N	Location	Pulling_C	No_of_Pul	Pulling_Req		
2	1	ma	kim	12-1-12-1	2011-01-C	1	Lee		
3									
4									
5									
6									

- Clicking "Export to Excel" will automatically be saved as "Pulling Record + Pulling Date.xls" in SBMS2.0/EXCEL folder.

Prescription Dept.	Prescription Dr.	Pulling Record	Total No. of Sam...
DP	kim	N	6
DR	Lee	Y	5


- When segment pulling is executed, "Pulling Record" values in Sample Databases change from "N" to "Y"

## 7) Export to Excel

A1																			
	A	B	C	D	E	F	G	H	I	J	K	L	M	N	O	P	Q	R	S
1	Sample_No	Patient_Name	Age	Sex	Registrat	Sample_Ty	Diagnosis	No_Remain	Location	Prescript	Prescript	Pulling_F	Total_No_	Total_Sam	Remark	Comment	Personal_ID		
2	6	sue	40	M	2011-01-CPH	DO	?	1-1-1-a11DA	kim	N	?	1ml					927291271928		
3	5	smun	29	M	2011-01-CPH	DO	?	1-1-1-a11DA	kim	N	?	1ml					548133291021		
4	4	stu	29	M	2011-01-CPH	DO	?	1-1-1-a11DA	park	N	?	1ml					548131164221		
5	3	master	29	M	2011-01-CPH	DO	?	1-1-1-a11DA	park	N	?	1ml					548133164971		
6	2	dai	13	M	2011-01-CRA	DA	6	3-1-1-231DP	kim	N	?	6	1ml	Test	Test		123452102893		
7	1	ma	12	F	2011-01-CRA	DA	1	12-1-12-1DR	Lee	Y	?	5	1ml	Test	Test		123456123456		
8																			
9																			
10																			
11																			
12																			
13																			
14																			
15																			
16																			
17																			
18																			
19																			
20																			

- Click "Export" from Main Menu to transfer the databases in the Archive to Microsoft Excel Spreadsheet
  - Confirm the data transferred to the Spreadsheet
- In the "Pulling Record" window, clicking "Export to Excel" will open the Microsoft Excel with the pulling records. In addition, there is also a function to export corresponding databases to Excel Spreadsheet in Main Frame.
  - Clicking "Export" will also automatically be saved as "Sample Information + Current Date/Time.xls" in SBMS2.0/EXCEL folder.

## 8) Configuration

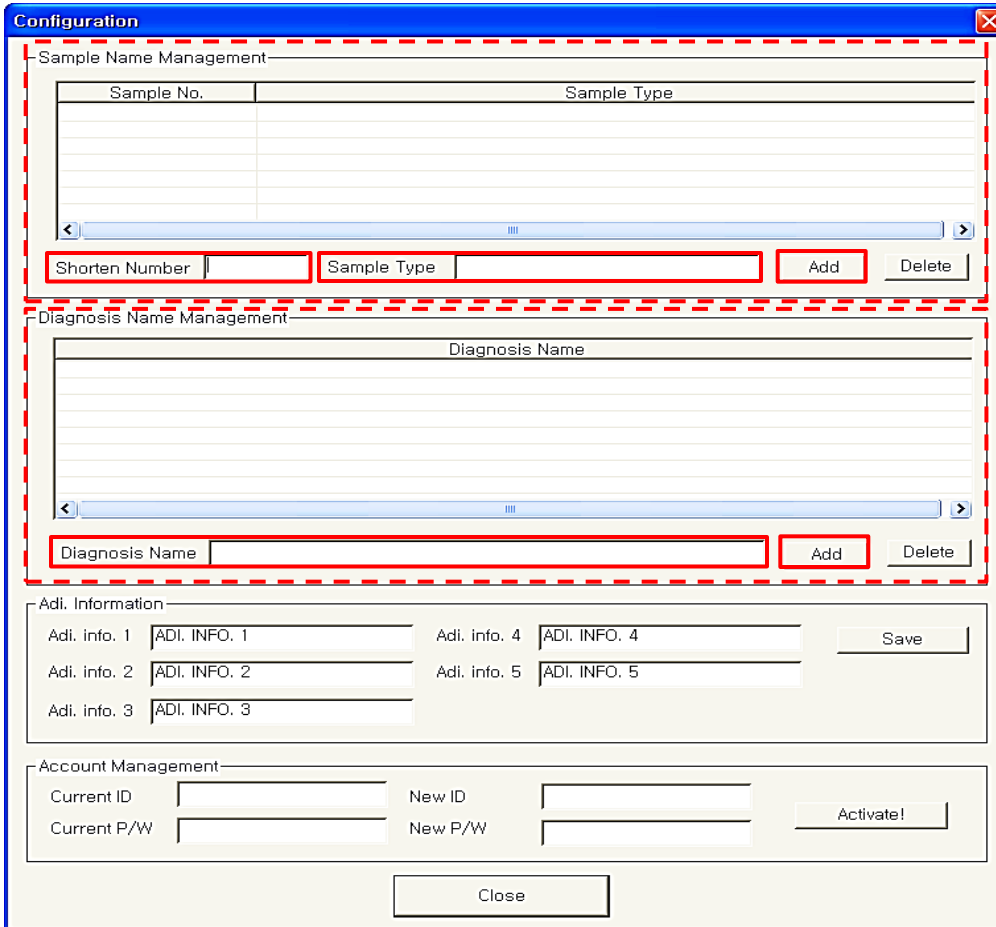

The screenshot shows a 'Configuration' window with the following sections:

- Sample Name Management:** A table with columns 'Sample No.' and 'Sample Type'. Below the table are input fields for 'Shorten Number' and 'Sample Type', and 'Add' and 'Delete' buttons.
- Diagnosis Name Management:** A table with a single column 'Diagnosis Name'. Below the table is an input field for 'Diagnosis Name' and 'Add' and 'Delete' buttons.
- Adi. Information:** Five input fields labeled 'Adi. info. 1' through 'Adi. info. 5', each containing the text 'ADI. INFO. X'. A 'Save' button is to the right.
- Account Management:** Input fields for 'Current ID', 'Current P/W', 'New ID', and 'New P/W'. An 'Activate!' button is to the right.

A 'Close' button is located at the bottom center of the window.


- ① Click "Config" from Main Menu to open "Configuration" window

### Assigning Sample Types

- ① Fill in "Sample Type" category with a title for the sample type
- ② Put a number to represent the sample type in "Shorten Number" category
- ③ Click "Add" to assign a sample type

### Assigning Diagnosis Names

- ① Fill in "Diagnosis Name" category with a title for the diagnosis
- ② Click "Add" to assign a diagnosis name

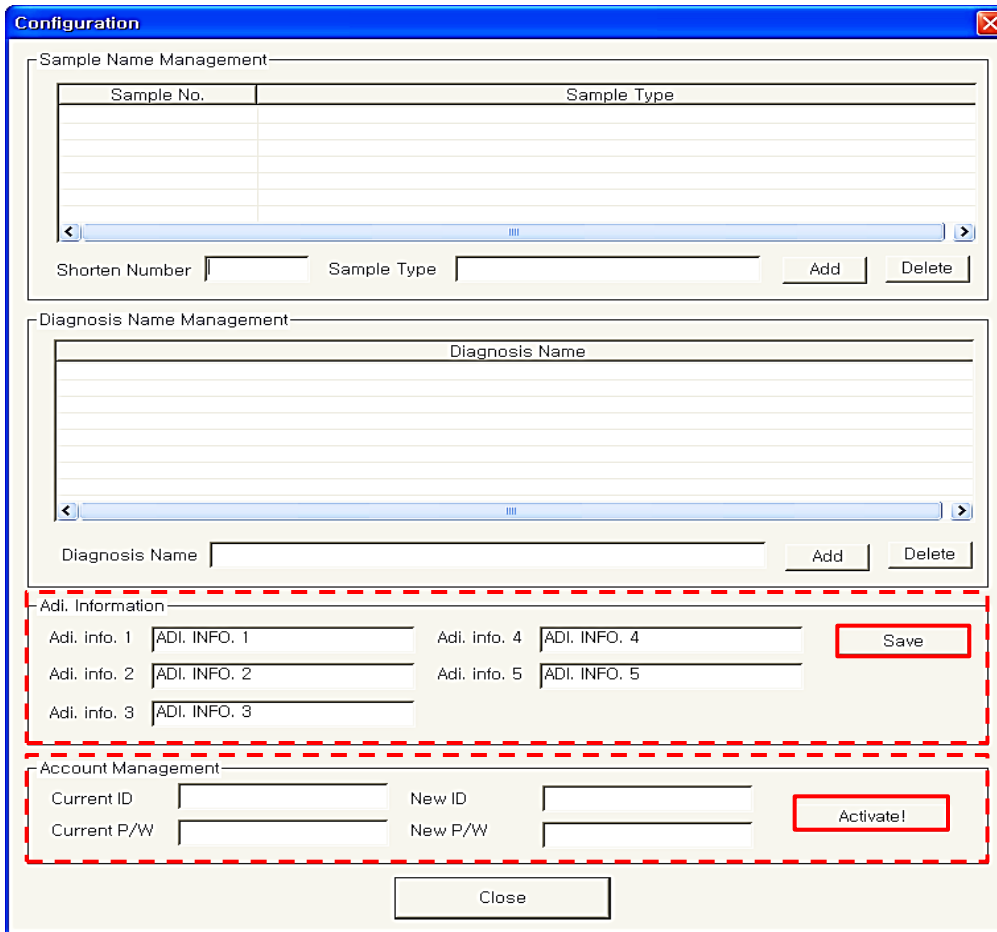


### Setup Additional Information


- ① Fill in any "Adi. info. #" under "Adi. Information" with a title for the additional information
- ② Click "Save" to assign an additional category


### Create or change current ID and Password

- ① Type current ID in "Current ID" and current Password in "Current P/W" under "Account Management"
- ② Fill in "New ID" and "New P/W"
- ③ Click "Activate!"
  - If there is no current ID or P/W, leave "Current ID" or "Current P/W" category blank.

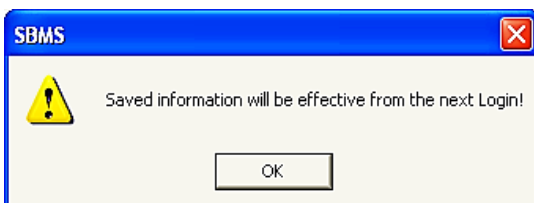

- New ID and Password will be effective on next login.
- For setting no ID or Password, leave "New ID" or "New P/W" category blank.

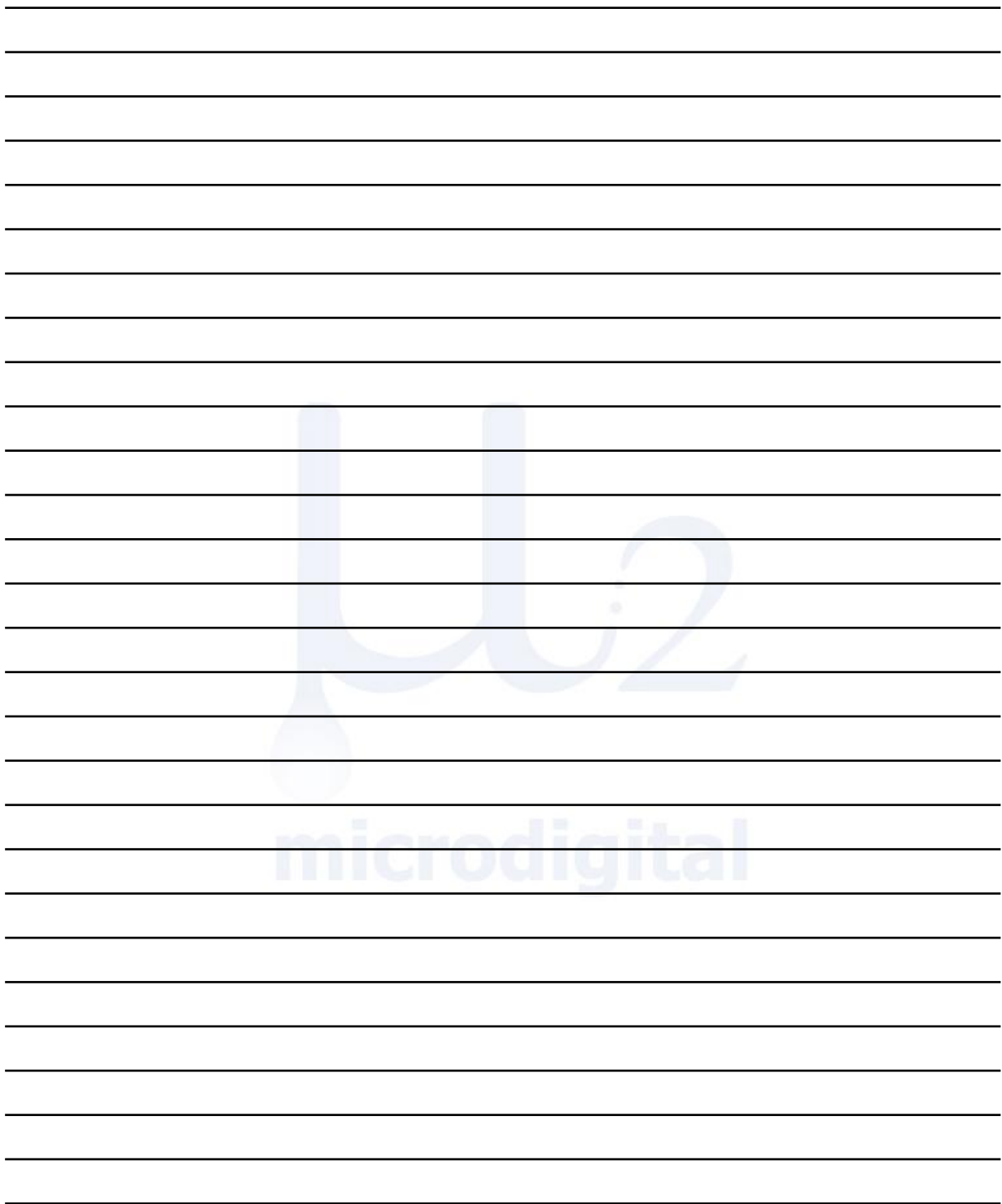

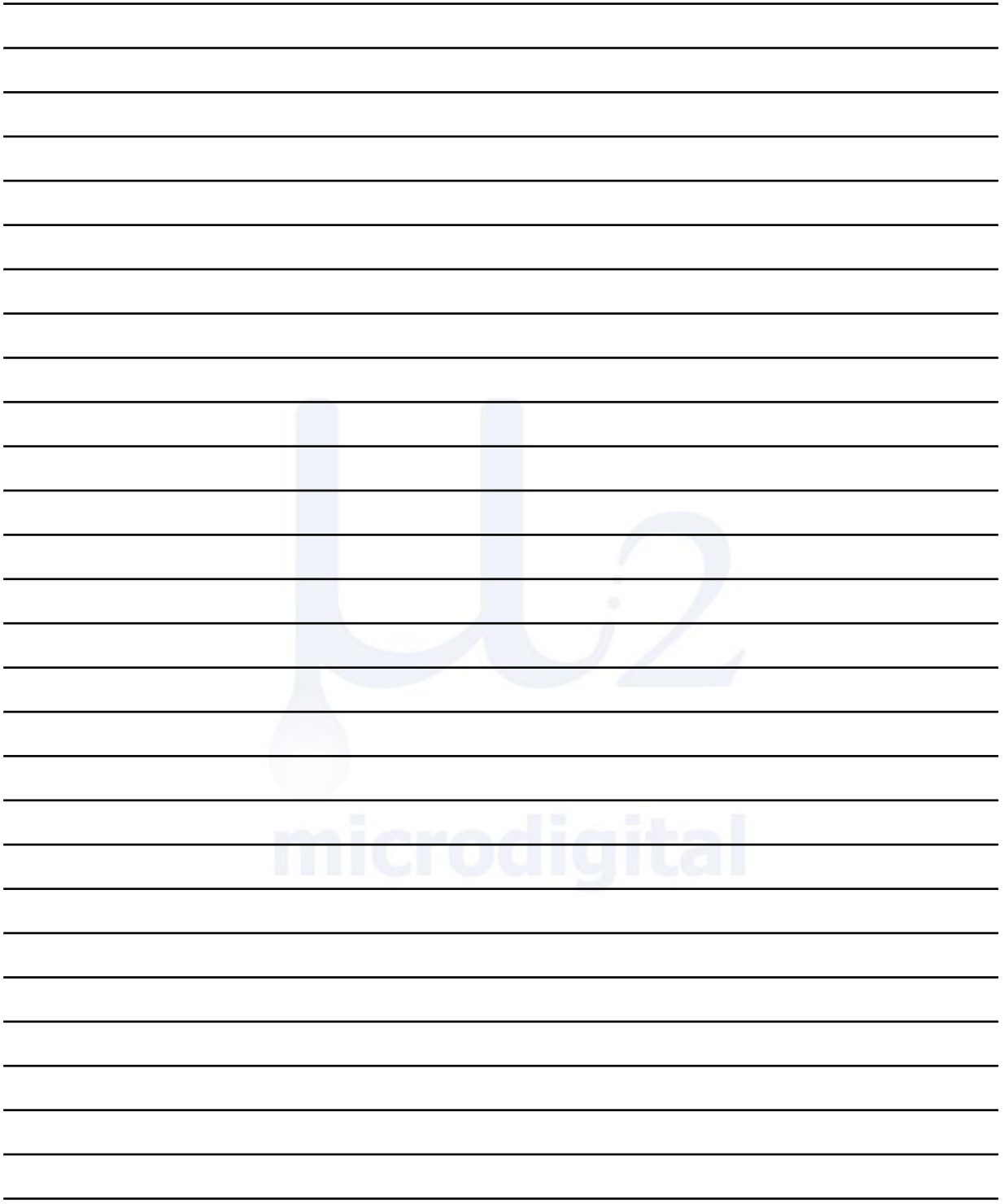

# $\mu$ 2

microdigital



**MicroDigital Co., Ltd.**  
#A-101, Korea Bio Park B/D, 694-1, SamPyung-Dong,  
BunDang-Gu, SungNam-Si, GyungGi-Do, 463-400, Korea  
Tel : +82-31-628-2200 / Fax : +82-31-628-2222  
[info@md-best.com](mailto:info@md-best.com) / [www.MD-Best.com](http://www.MD-Best.com)