# RayBio<sup>®</sup> Human Cytokine Antibody Array G Series 4000

### **Patent Pending Technology**

#### User Manual (Revised February 18, 2009)

Combination of Human Cytokine Antibody Array G Series 6, 7, 8, 9 & 10
Cat# AAH-CYT-G4000-4

RayBio<sup>®</sup> Human Cytokine Antibody Array G Series 6 (Cat# AAH-CYT-G6-4) RayBio<sup>®</sup> Human Cytokine Antibody Array G Series 7 (Cat# AAH-CYT-G7-4) RayBio<sup>®</sup> Human Cytokine Antibody Array G Series 8 (Cat# AAH-CYT-G8-4)

RayBio<sup>®</sup> Human Cytokine Antibody Array G Series 9(Cat# AAH-CYT-G9-4) RayBio<sup>®</sup> Human Cytokine Antibody Array G Series 10 (Cat# AAH-CYT-G10-4

Combination of Human Cytokine Antibody Array G Series 6, 7, 8, 9 & 10 Cat# AAH-CYT-G4000-8

RayBio<sup>®</sup> Human Cytokine Antibody Array G Series 6 (Cat# AAH-CYT-G6-8) RayBio<sup>®</sup> Human Cytokine Antibody Array G Series 7 (Cat# AAH-CYT-G7-8) RayBio<sup>®</sup> Human Cytokine Antibody Array G Series 8 (Cat# AAH-CYT-G8-8) RayBio<sup>®</sup> Human Cytokine Antibody Array G Series 9(Cat# AAH-CYT-G9-8) RayBio<sup>®</sup> Human Cytokine Antibody Array G Series 10 (Cat# AAH-CYT-G10-8)

Please read the manual carefully before you start your experiment



We Provide You With Excellent Protein Array Systems And Service

Tel:(Toll Free) 1-888-494-8555 or 770-729-2992; Fax: 1-888-547-0580; Website:www.raybiotech.com Email: info@raybiotech.com



RayBio® Human Cytokine Antibody Array G Series 4000 Protocol

## **TABLE OF CONTENTS**

I.	Introduction	2
	How It Works	5
II.	Materials Provided	6
	Additional Materials Required	6
III.	Overview and General Considerations	7
	A. Preparation of Samples	7
	B. Handling Glass Chips	8
	C. Incubation.	8
IV.	Protocol	9
	A. Blocking and Incubation	9
	B. Detection.	10
V.	Interpretation of Results	13
VI.	Troubleshooting Guide	21
VII.	Selected References Using RayBiotech Products	22

Cytokine Antibody Arrays are RayBiotech patent-pending technology.

RayBio<sup>®</sup> is the trademark of RayBiotech, Inc.

#### I. Introduction

All cell functions, including cell proliferation, cell death and differentiation, as well as maintenance of health status and development of disease, are controlled by many genes and signaling pathways. New techniques such as cDNA microarrays have enabled us to analyze the global gene expression <sup>1-3</sup>. However, almost all cell functions are executed by proteins, which cannot be studied by DNA and RNA alone. Experimental analysis clearly shows a disparity between the relative expression levels of mRNA and their corresponding proteins <sup>4</sup>. Therefore, it is critical to analyze the protein profile. Currently, two-dimensional polyacrylamide SDS page coupled with mass spectrometry is the mainstream approach to analyzing multiple protein expression levels <sup>5,6</sup>. However, the requirement of sophisticated devices and the lack of quantitative measurements greatly limit its broad application. Thus, no simple, cost effective, and rapid method of analysis of multiple protein expression levels has been available to researchers until now.

Our RayBio<sup>®</sup> Human Cytokine Antibody Array is the first commercially available protein array system <sup>7-11</sup>. By using the RayBiotech system, scientists can rapidly and accurately identify the expression profiles of multiple cytokines in several hours inexpensively.

The RayBiotech kit (G series) is a glass chip format. The kit provides a highly sensitive approach to simultaneously detect multiple cytokine expression levels from cell culture supernatant, patient's serum, tissue lysate and other sources. The arrays are manufactured using non-contact arrayer. The experimental procedure is simple and can be performed in any laboratory. The signals from G series arrays are detected using a laser scanner.

The RayBio<sup>®</sup> Human Cytokine Antibody Array G series 4000 is the only product available in the market that can detect 274 human cytokines in single experiment.

Pathway-specific array systems allow investigators to focus on the specific problem and are becoming an increasingly powerful tool in cDNA microarray system. RayBiotech's first protein array system, known as

RayBio<sup>®</sup> Human Cytokine Antibody Array, is particularly useful compared with the human cytokine cDNA microarray system. Besides the ability to detect protein expression, RayBiotech's system is a more accurate reflection of active cytokine levels because it only detects secreted cytokines, and no amplification step is needed. Furthermore, it is much simpler, faster, environmentally friendlier, and more sensitive.


Simultaneous detection of multiple cytokines undoubtedly provides a powerful tool to study cytokines. Cytokines play an important role in innate immunity, apoptosis, angiogenesis, cell growth and differentiation <sup>12</sup>. Cytokines are involved in most disease processes, including cancer and cardiac diseases. The interaction between cytokines and the cellular immune system is a dynamic process. The interactions of positive and negative stimuli, and positive as well as negative regulatory loops are complex and often involve multiple cytokines.

#### **Reference List**

- 1. LPS induces the interaction of a transcription factor, LPS-induced TNF-a factor, and STAT6(B) with effects on multiple cytokines. Xiaoren Tang, Deborah Levy Marciano, Susan E. Leeman and Salomon Amar. **PNAS**. April 5, 2005 vol. 102 no. 14 5132-5137
- 2. HIV-1-mediated apoptosis of neuronal cells: Proximal molecular mechanisms of HIV-1-induced encephalopathy. Yan Xu, Joseph Kulkoshy, Roger j. Pomerantz. **PNAS**. 2004 May 4, 2004 Vol. 101 No. 18.
- 3. Synergistic increases in intracellular Ca(2+), and the release of MCP-1, RANTES, and IL-6 by astrocytes treated with opiates and HIV-1 Tat. **GLIA**. 2005 Apr 15;50(2):91-106.
- 4. Bone Marrow Stroma Influences Transforming Growth Factor-β Production in Breast Cancer Cells to Regulate c-myc Activation of the Preprotachykinin-I Gene in Breast Cancer Cells. Hyun S. Oh, Anabella Moharita... Pranela Rameshwar. **CANCER RESEARCH**. 64, 6327-6336, September 1, 2004

- 5. Recombinant Herpes Simplex Virus Type 1 (HSV-1) Codelivering Interleukin-12p35 as a Molecular Adjuvant Enhances the Protective Immune Response against Ocular HSV-1 Challenge. **J Virology**. Mar. 2005 Vol. 79, No. 6.
- 6. Dysregulated Inflammatory Response to *Candida albicans* in a C5-Deficient Mouse Strain. Alaka Mullick, Miria Elias, Serge Picard, Philippe Gros. **Infection and Immunity**, Oct. 2004, p. 5868-5876, DOI: 10.1128/IAI.72.10.5868-5876.2004
- 7. Leukotriene B<sub>4</sub> Strongly Increases Monocyte Chemoattractant Protein-1 in Human Monocytes Li Huang, Annie Zhao, Frederick Wong, Julia M. Ayala, Jisong Cui **Arteriosclerosis, Thrombosis, and Vascular Biology**. 2004;24:1783-1788
- 8. Human CD1d-unrestricted NKT cells release chemokines upon Fas engagement. Martin Giroux and François Denis. Yan Xu, Joseph Kulkoshy, Roger j. Pomerantz. **Blood**. prepublished online September 2, 2004; DOI 10.1182/blood-2004-04-1537
- 9. Monitoring the response of orthotopic bladder tumors to granulocyte macrophage colony-stimulating factor therapy using the prostate-specific antigen gene as a reporter. Wu Q, Esuvaranathan K, Mahendran R. **Clinical Cancer Research**. 2004 Oct 15; 10(20):6977-84.
- 10. Neuroglial activation and neuroinflammation in the brain of patients with autism (p NA). Diana L. Vargas, Caterina Nascimbene, Chitra Krishnan, Andrew W. Zimmerman, Carlos A. Pardo. **Annals of Neurology.** 2005 Jan 1; DOI: 10.1002/ana.20315
- 11. Cytokine profiling of macrophages exposed to Porphyromonas gingivalis, its LPS or its FimA. Qingde Zhou, Tesfahun Desta, Dana T. Graves and Salomon Amar. **Infection and Immunity** (IAI). 2005 Feb;73(2):935-43.
- 12. Veto-like activity of mesenchymal stem cells: functional discrimination between cellular responses to alloantigens and recall antigens. Rameshwar P. **Journal of Immunology**. 2003 Oct 1;171(7):3426-34.
- 13. The promise of cytokine antibody arrays in drug discovery process. Huang RP, Yang W, Yang Y, Flowers L, et al. **Expert Opinion on Drug Discovery**. (2005) 9:601-615.

### Here's how it works

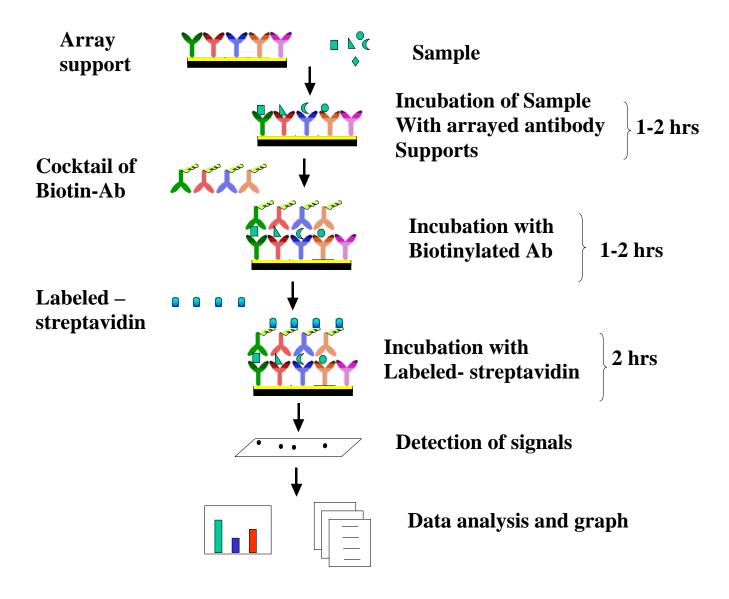

#### **II. Materials Provided**


Upon receipt, all components of the RayBio<sup>®</sup> Human Cytokine Antibody Array kit should be stored at -20°C. At -20°C the kit will retain complete activity for up to 6 months. Once thawed, the glass chips, Fluorescent dyestreptavidin, Internal Control and 2X Blocking Buffer should be kept at –20°C and all other component should be stored at 4°C. Use within three months after reagents have been thawed. Please use within six months of purchase.

- RayBio<sup>®</sup> Human Cytokine Antibody Microarray slides (5 slides with 4 or 8 subarrays each: Human 6, 7, 8, 9, and 10)
- Biotin-Conjugated Anti-Cytokines (1 tube/4 subarrays, 1/2 tubes each Human 6, 7, 8, 9, and 10)
- 1,500X Fluorescent Dye-conjugated Streptavidin (Cy3 equivalent, 1 tube per 20 subarrays, 1 or 2 tubes)
- 2X Blocking Buffer (10 ml)
- 20X Wash Buffer I (30 ml X 2)
- 20X Wash Buffer II (30 ml X 2)
- Internal control (powder, 1 tube)
- 2X Cell Lysis Buffer (10 ml)
- RayBio<sup>®</sup> G series antibody array accessory (including slide incubation chamber, Gasket, Protective cover, Snap-on sides and adhesive film)
- 30 ml tube
- Manual


### **Additional Materials Required**

- Orbital shaker
- Laser scanner for fluorescence detection
- Aluminum foil
- Distilled water
- Plastic box

## Layout of G series 4000

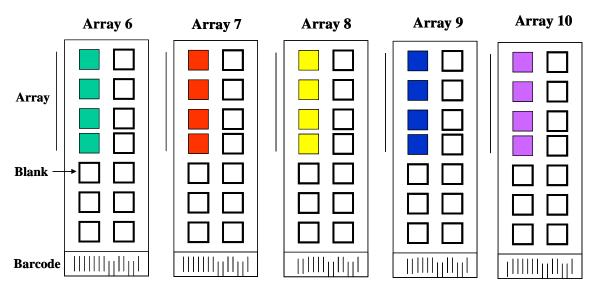

4 arrays in one glass chip

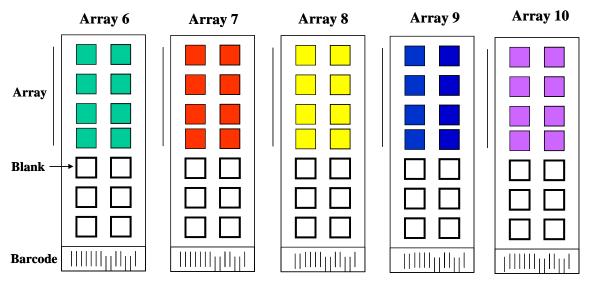

8 arrays in one glass chip

## III. Overview and General Considerations

## A. Preparation of Samples

- Use serum-free conditioned media if possible.
- If serum-containing conditioned media is required, use uncultured media as a negative control sample, since many types of sera contain cytokines.

- For cell lysates and tissue lysates, we recommend using RayBio® Cell Lysis Buffer to extract proteins from cell or tissue (e.g. using homogenizer). Dilute 2X RayBio® Cell Lysis Buffer with H<sub>2</sub>O (we recommend adding proteinase inhibitors to Cell Lysis Buffer before use). After extraction, spin the sample down and save the supernatant for your experiment. Determine protein concentration.
- We recommend using:
  - o 50–100 μl of Conditioned media (undiluted), or
  - o 50–100 μl of 2-fold to 5-fold diluted sera or plasma, or
  - o 10–200 μg of total protein for cell lysates and tissue lysates.

If you experience high background, you may further dilute your sample.

#### B. Handling glass chips

- The microarray slides are sensitive, do not touch the surface. Hold the slides by the edges only.
- Handle all buffers and slides with powder-free gloves.
- Avoid breaking glass slide.
- Handle glass chip in clean environment.

### C. Incubation


- Completely cover array area with sample or buffer during incubation, and cover the incubation chamber with adhesive film or plastic sheet protector to avoid drying.
- Avoid foaming during incubation steps.
- Perform all incubation and wash steps under gentle rotation.
- Cover the incubation chamber with adhesive film during incubation, particularly when incubation is more than 2 hours or 50 µl of sample or reagent is used.
- Avoid cross-contamination from overflowing solution to neighboring wells.
- Several incubation steps such as step 3 (blocking), step 4 (sample incubation), step 9 (biotin-Ab incubation) or step 12 (Fluorescent dyestreptavidin incubation) may be done at 4°C for overnight. Please make sure to cover the incubation chamber tightly to prevent evaporation.


#### IV. Protocol

### A. Blocking and Incubation

- 1. Take the glass chip out from the box. Let air dry for 2 hours.
- 2. Assemble the glass chip into incubation chamber and incubation frame as shown below. (Note: if you slide has be assembled, you can go to step 3 directly).



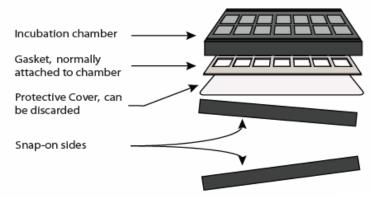

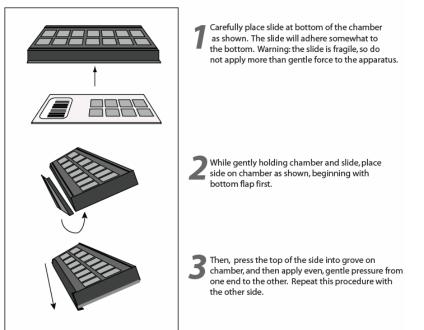

3. Add 100 µl 1 X Blocking Buffer into each well and incubate at room temperature for 30 min to block slides. Dilute 2X Blocking Buffer with H<sub>2</sub>O. Make sure no bubbles are in the well.

*Note: only add reagents to wells printed with antibodies.* 

4. Decant Blocking Buffer from each well, Add 50 to 100 μl of each sample to subarray 6, subarray 7, subarray 8, subarray 9, and subarray 10. Incubate arrays with sample at room temperature for 1 to 2 hours. Dilute sample using 1X Blocking Buffer if necessary. We strongly recommend including Internal Control (IC) in your assay. Add 100 μl of Blocking Buffer to IC tube, mix well and transfer 1 μl of IC to each well (50 to 100 μl of sample).

Note: when you transfer IC, use 0.1 µl to 2 µl pipettor.

*Note: Incubation may be done at*  $4^{\circ}C$  *for overnight.* 

Note: We recommend using 50 to 100 µl of conditioned media or 50 to 100 µl of 2-5 fold diluted serum or plasma or 10-200 µg of protein for cell lysates and tissue lysates. Dilute the lysate at least 10 fold with 1 X blocking buffer to make a total volume of 50 to 100 µl. Make sure there are no bubbles in the wells.

Note: The amount of sample used depends on the abundance of cytokines. More concentrated sample can be used if signals are too weak. If signals are too strong, the sample can be diluted further.

5. Decant the samples from each well, and wash 3 times with 150 μl of 1X Wash Buffer I at room temperature with gentle shaking. 2 min per wash. Dilute 20X Wash Buffer I with H<sub>2</sub>O. Completely remove wash buffer I in each wash step.

*Note: avoid solution flowing into neighboring wells.* 

- 6. Put the glass chip with frame into a box with 1 X Wash Buffer I (cover the whole glass slide and frame with Wash Buffer I), and wash 2 times at room temperature with gentle shaking for 10 min per each wash.
- 7. Decant the 1X Wash Buffer I from each well, Put the glass chip with frame into the box with 1X Wash Buffer II (cover the whole glass slide and frame with Wash Buffer II), and wash 2 times at room temperature

- with gentle shaking for 5 min per each wash. Remove all of Wash Buffer II in the well. Dilute 20X Wash Buffer II with H<sub>2</sub>O.
- 8. Prepare working solution for biotin-conjugated antibodies. After brief spin to force antibody aliquot to bottom of tube,
- A. Add 300 µl of 1X blocking buffer to the Biotin-Conjugated Antibody 6 tube. Mix gently.
- B. Add 300 μl of 1X blocking buffer to the Biotin-Conjugated Antibody 7 tube. Mix gently.
- c. Add 300 µl of 1X blocking buffer to the Biotin-Conjugated Antibody 8 tube. Mix gently.
- D. Add 300 µl of 1X blocking buffer to the Biotin-Conjugated Antibody 9 tube. Mix gently.
- E. Add 300 μl of 1X blocking buffer to the Biotin-Conjugated Antibody 10 tube. Mix gently.

*Note: the diluted biotin-conjugated antibodies can be stored at*  $4^{0}$ *C for* 2-3 *days.* 

9. Add 70 μl of diluted biotin-conjugated antibodies to each corresponding well (70 μl of diluted biotin-conjugated antibodies 6 to each subarray 6, 70 μl of diluted biotin-conjugated antibodies 7 to each subarray 7, 70 μl of diluted biotin-conjugated antibodies 8 to each subarray 8), 70 μl of diluted biotin-conjugated antibodies 9 to each subarray 9) and 70 μl of diluted biotin-conjugated antibodies 10 to each subarray 10). Incubate at room temperature for 2 hours.

*Note: incubation may be done at*  $4^{\circ}$ *C for overnight.* 

10. Wash as directed in steps 5 and then wash 3 times with 150 μl of 1X Wash Buffer II at room temperature with shaking. 2 min per wash. Completely remove wash buffer II in each wash step.

- 11. Add 70 µl of 1,500 fold diluted Fluorescent dye-conjugated streptavidin (after brief spinning, add 1.5 ml of Blocking Buffer to Fluorescent dye-conjugated streptavidin tube) to each subarray. Cover the incubation chamber with Adhesive film. Cover the plate with aluminum foil to avoid exposure to light or incubate in dark room.
- 12. Incubate at room temperature for 1 to 2 hours.

*Note: incubation may be done at 4°C for overnight.* 

13. Wash with Wash Buffer I **twice** as directed in steps 5.

#### **B.** Fluorescence Detection

- 1. Decant excess Wash Buffer from wells.
- 2. Disassemble the slide out of the incubation frame and chamber.
- 3. Place the whole slide in a 30 ml centrifuge tube provided, add enough Wash Buffer I (about 20 ml) to cover the whole slide and gently shake at room temperature for 10 minutes. Decant Wash Buffer I. Repeat Wash Buffer I once. Wash with Wash Buffer II (about 20 ml) with gentle shake at room temperature for 10 minutes. Or wash using slide chamber. Rinse the slide with distilled H<sub>2</sub>O.
- 4. Remove water droplets by centrifuge at 1,000 rpm for 3 minutes and then let slide dry completely in air at least 20 minutes (protect from light). Make sure the slides are absolutely dry before the scanning procedure.
- 5. Image the signals using laser scanner such as Axon GenePix using cy3 or "green" channel (Excitation frequency 532 nm).

Note: we recommend scanning slides right after experiment. You also can store the slide at  $-20^{\circ}$ C in dark for several days. If you do not have a laser scanner, RayBiotech can provide service for you. Just simply send your slide to us and we will take care of it.


### V. Interpretation of Results:

The following figure shows RayBio<sup>®</sup> Human Cytokine Antibody Array G series 2000 probed with different cell culture supernatant. The images were captured using laser scanner. The biotin-conjugated protein produces positive signals, which can be used to identify the orientation and to compare the relative expression levels among the different wells. The internal control (IC) can also be used to normalize the signal intensities among array membranes in different experiments.

The signal intensities obtained from laser scanner can simply be imported into our analysis tool. The analysis tool will help you:

- Locate your signal intensities to antibody array map
- Link the protein to website for more detailed information on the particular protein
- Protein list sorting
- Average signal intensities
- Subtract background
- Normalize the data from different samples
- Obtain protein level comparison charts among different samples

This analysis tool is very simple and affordable, which will not only assist in compiling and organizing your data, but also reduces your calculations to a "copy and paste" step.

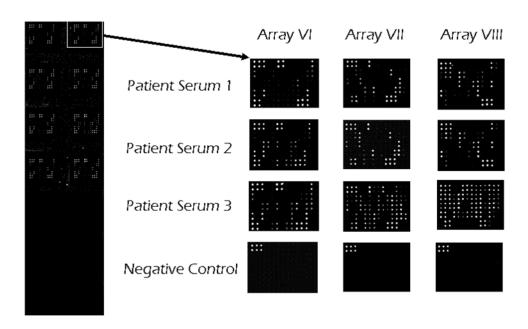

If you do not use our **RayBio**<sup>®</sup> **Analysis Tool,** you can locate the cytokines by referring to RayBio<sup>®</sup> Human cytokine Antibody Array G series 2000. Please keep in mind that G series 2000 consists of three individual arrays; human cytokine antibody array 6, human cytokine antibody array 7 and human cytokine antibody array 8. Refer to corresponding map.

#### Normalization and comparison

For biomarker discovery or for analysis of large number of arrays, great attention must be paid to the normalization. Our antibody array design includes several controls for normalization and comparison of arrays performing in different membranes and different experiments (for more information please read the reference 13 from the introduction).

**Positive control.** Positive control is biotinylated protein. It can be used to normalize the streptavidin incubation step. If the positive signals from different array membranes are similar, positive control is a simple and effective way for normalization.

*Internal control*. RayBio<sup>®</sup> antibody arrays also include spiking-in protein serving as internal control (IC). The spiking-in proteins do not have cross-reactivity with protein in the array. It can be used to normalize the entire process.

*Negative control*. Negative control is BSA. Normally, it should only give a background reading.

#### **Data Extraction Tips:**

- Ignore any comet tails
- Define the area for signal capture for all spots as 110-120 micron diameter, using the same area for every spot.
- Use median signal value, not the total or the mean
- Use local background correction (also median value).
- Exclude obvious outlier data in its calculations.

Using these guidelines, along with using PMT, brightness and contrast settings that reduce the background as much as possible, we get very good results with interassay and intraassay CV <20%, even with some imperfections in the antibody spots.

Threshold of significant difference in expression: Any  $\geq 1.5$ -fold increase or  $\leq 0.65$ -fold decrease in signal intensity for a single analyte between samples, provided that both signals are well above background (Mean background + 2 standard deviations, accuracy  $\approx 95\%$ ).

#### Array Map Notes:

- o IL-12 reacts with both IL-12p40 and IL-12p70.
- o IL-12p70 only recognizes IL-12p70.
- o GRO reacts with CXCL1, CXCL2 and CXCL3 (GRO alpha, beta, and gamma, respectively).
- o GRO alpha reacts only with CXCL1.
- VEGF reacts with VEGF-165 and VEGF-121
- o TGF-beta 1 reacts only with active form of TGF-beta 1
- o All MMP antibody pairs detect both pro and mature forms

Abbreviations: IP-10, Interferon-inducible protein-10; LAP, latency associated peptide (TGF-β1); LH, leutinizing hormone; LIF, leukocyte inhibitory factor. MMP, Matrix Metalloproteinase; PF, plasma factor 4; Pos, positive control; Neg, negative control. All other are used standard abbreviations.

## RayBio® Human Cytokine Antibody Array G series 6 (60)

	Α	В	С	D	E	F	G	Н	ı	J	К	L	М	N
1	POS 1	POS 2	POS 3	NEG	NEG	Angiogenin	BDNF	BLC	BMP-4	BMP-6	CK b 8-1	CNTF	EGF	Eotaxin
2	POS 1	POS 2	POS 3	NEG	NEG	Angiogenin	BDNF	BLC	BMP-4	BMP-6	CK b 8-1	CNTF	EGF	Eotaxin
3	Eotaxin-2	Eotaxin-3	FGF-6	FGF-7	Fit-3 Ligand	Fractalkine	GCP-2	GDNF	GM-CSF	I-309	IFN- gamma	IGFBP-1	IGFBP-2	IGFBP-4
4	Eotaxin-2	Eotaxin-3	FGF-6	FGF-7	Fit-3 Ligand	Fractalkine	GCP-2	GDNF	GM-CSF	1-309	IFN- gamma	IGFBP-1	IGFBP-2	IGFBP-4
5	IGF-I	IL-10	IL-13	IL-15	IL-16	IL-1alpha	IL-1beta	IL-1ra	IL-2	IL-3	IL-4	IL-5	IL-6	IL-7
6	IGF-I	IL-10	IL-13	IL-15	IL-16	IL-1alpha	IL-1beta	IL-1ra	IL-2	IL-3	IL-4	IL-5	IL-6	IL-7
7	Leptin	LIGHT	MCP-1	MCP-2	MCP-3	MCP-4	M-CSF	MDC	MIG	MIP-1 delta	MIP-3 alpha	NAP-2	NT-3	PARC
8	Leptin	LIGHT	MCP-1	MCP-2	MCP-3	MCP-4	M-CSF	MDC	MIG	MIP-1 delta	MIP-3 alpha	NAP-2	NT-3	PARC
9	PDGF-BB	RANTES	SCF	SDF-1	TARC	TGF-beta1	TGF-beta 3	TNF-alpha	TNF-beta	IC1	IC2	IC3	NEG	NEG
10	PDGF-BB	RANTES	SCF	SDF-1	TARC	TGF-beta1	TGF-beta 3	TNF-alpha	TNF-beta	IC1	IC2	IC3	NEG	NEG

# RayBio® Human Cytokine Antibody Array G series 7 (60)

	Α	В	С	D	E	F	G	Н		J	К	L	М	N
1	POS1	POS 2	POS 3	NEG	NEG	Acrp30	AgRP	Angio- poietin-2	Amphi- regulin	AxI	bFGF	b-NGF	втс	CCL-28
2	POS1	POS 2	POS 3	NEG	NEG	Acrp30	AgRP	Angio- poietin-2	Amphi- regulin	AxI	bFGF	b-NGF	втс	CCL-28
3	CTACK	Dtk	EGF-R	ENA-78	Fas/ TNFRSF6	FGF-4	FGF-9	GCSF	GITR-Ligand	GITR	GRO	GRO- alpha	HCC-4	HGF
4	CTACK	Dtk	EGF-R	ENA-78	Fas/ TNFRSF6	FGF-4	FGF-9	GCSF	GITR-Ligand	GITR	GRO	GRO- alpha	HCC-4	HGF
5	ICAM-1	ICAM-3	IGFBP-3	IGFBP-6	IGF-I SR	IL-1 R4/ST2	IL-1 RI	IL-11	IL-12 p40	IL-12 p70	IL-17	IL-2 R alpha	IL-6 R	IL-8
6	ICAM-1	ICAM-3	IGFBP-3	IGFBP-6	IGF-I SR	IL-1 R4/ST2	IL-1 RI	IL-11	IL-12 p40	IL-12 p70	IL-17	IL-2 R alpha	IL-6 R	IL-8
7	I-TAC	Lympho- tactin	MIF	MIP-1 alpha	MIP-1 beta	MIP-3 beta	MSP-alpha	NT-4	Osteo- protegerin	Oncostatin M	PIGF	sgp130	sTNF RII	sTNF-RI
8	I-TAC	Lympho- tactin	MIF	MIP-1 alpha	MIP- 1beta	MIP-3 beta	MSP-alpha	NT-4	Osteo- protegerin	Oncostatin M	PIGF	sgp130	sTNF RII	sTNF-RI
9	TECK	TIMP-1	TIMP-2	Thrombo- poietin	TRAIL R3	TRAIL R4	uPAR	VEGF	VEGF-D	IC 1	IC 2	IC 3	NEG	NEG
10	TECK	TIMP-1	TIMP-2	Thrombo- poietin	TRAIL R3	TRAIL R4	uPAR	VEGF	VEGF-D	IC 1	IC 2	IC 3	NEG	NEG

# RayBio® Human Cytokine Antibody Array G series 8 (54)

	Α	В	С	D	Ε	F	G	н	I	J	K	L	М	N
1	POS 1	POS 2	POS 3	NEG	NEG	Activin A	ALCAM	B7-1 (CD80)	BMP-5	BMP-7	Cardio- trophin-1	CD14	CXCL- 16	DR6 (TNFRSF21)
2	POS 1	POS 2	POS 3	NEG	NEG	Activin A	ALCAM	B7-1 (CD80)	BMP-5	BMP-7	Cardio- trophin-1	CD14	CXCL- 16	DR6 (TNFRSF21)
3	Endoglin	ErbB3	E-Selectin	Fas Ligand	ICAM-2	IGF-II	IL-1 R II	IL-10 R beta	IL-13 R alpha 2	IL-18 BP alpha	IL-18 R beta	MMP-3	IL-2 R beta	IL-2 R gamma
4	Endoglin	ErbB3	E-Selectin	Fas Ligand	ICAM-2	IGF-II	IL-1 R II	IL-10 R beta	IL-13 R alpha 2	IL-18 BP alpha	IL-18 R beta	MMP-3	IL-2 R beta	IL-2 R gamma
5	IL-21R	IL-5 R alpha	IL-9	IP-10	LAP	Leptin R	LIF	L-Selectin	M-CSF R	MMP-1	MMP-13	MMP-9	MPIF-1	NGF R
6	IL-21R	IL-5 R alpha	IL-9	IP-10	LAP	Leptin R	LIF	L-Selectin	M-CSF R	MMP-1	MMP-13	MMP-9	MPIF-1	NGF R
7	PDGF AA	PDGF-AB	PDGF R alpha	PDGF R beta	PECAM-1	Prolactin	SCF R	SDF-1 beta	Siglec-5	TGF-alpha	TGF-beta 2	Tie-1	Tie-2	TIMP-4
8	PDGF AA	PDGF-AB	PDGF R alpha	PDGF R beta	PECAM-1	Prolactin	SCF R	SDF-1 beta	Siglec-5	TGF-alpha	TGF-beta 2	Tie-1	Tie-2	TIMP-4
9	VE- Cadherin	VEGF R2	VEGF R3	Neg	Neg	Neg	Neg	Neg	Neg	IC1	IC2	IC3	NEG	NEG
10	VE- Cadherin	VEGF R2	VEGF R3	Neg	Neg	Neg	Neg	Neg	Neg	IC1	IC2	IC3	NEG	NEG

# RayBio® Human Cytokine Antibody Array G series 9 (51)

	а	b	С	d	е	f	g	h	i	j	k	ĺ
1	POS1	POS2	POS3	NEG	NEG	Adiposin	BCAM	CD30	CD40	Fcr RIIB/C	Ferritin	FLRG
2	POS1	POS2	POS3	NEG	NEG	Adiposin	BCAM	CD30	CD40	Fcr RIIB/C	Ferritin	FLRG
3	Follistatin	Furin	Galectin-7	GDF-15	Growth Hormon	IL-10 R alpha	IL-22	IL-28A	IL29	IL-31	Insulin	Luteinizing hormone
4	Follistatin	Furin	Galectin-7	GDF-15	Growth Hormon	IL-10 R alpha	IL-22	IL-28A	IL29	IL-31	Insulin	Luteinizing hormone
5	LIMPII	LYVE-1	Marapsin	MICA	MICB	MMP-2	MMP-7	MMP-8	MMP-10	NCAM-1	Nidogen-1	NrCAM
6	LIMPII	LYVE-1	Marapsin	MICA	MICB	MMP-2	MMP-7	MMP-8	MMP-10	NCAM-1	Nidogen-1	NrCAM
7	NRG1-beta1	Osteopontin	PAI-I	Platelet factor 4	PSA-total	RAGE	RANK	Resistin	SAA	Siglec-9	TACE	TIM-1
8	NRG1-beta1	Osteopontin	PAI-I	Platelet factor 4	PSA-total	RAGE	RANK	Resistin	SAA	Siglec-9	TACE	TIM-1
9	TRAIL R2	Trappin-2	TREM-1	TSH	TSLP	VCAM-1	VEGF-C	XEDAR	NEG	IC 1	IC 2	IC 3
10	TRAIL R2	Trappin-2	TREM-1	TSH	TSLP	VCAM-1	VEGF-C	XEDAR	NEG	IC 1	IC 2	IC3

## RayBio® Human Cytokine Antibody Array G series 10 (49)

	а	b	С	d	е	f	g	h	i	j	k	ı
1	POS1	POS2	POS3	NEG	NEG	4-1BB	ACE-2	Alpha-fetoprotein	Angiopoietin-1	Angiostatin	ANGPTL4	Bate2 M
2	POS1	POS2	POS3	NEG	NEG	4-1BB	ACE-2	Alpha-fetoprotein	Angiopoietin-1	Angiostatin	ANGPTL4	Bate2 M
3	BCMA	beta IG-H3	CA125	CA15-3	CA19-9	Carbonic Anhydrase IX	Cathepsin S	CCL14a	CCL21	CD23	CD40 Ligand	CEA
4	BCMA	beta IG-H3	CA125	CA15-3	CA19-9	Carbonic Anhydrase IX	Cathepsin S	CCL14a	CCL21	CD23	CD40 Ligand	CEA
5	CEACAM-1	Cripto-1	CRP	DAN	Decorin	DKK-1	DKK-3	DKK-4	DPPIV	E-Cadherin	EDA-A2	EG-VEGF
6	CEACAM-1	Cripto-1	CRP	DAN	Decorin	DKK-1	DKK-3	DKK-4	DPPIV	E-Cadherin	EDA-A2	EG-VEGF
7	EpCAM	ErbB2	Erythropoietin R	FSH	HB-EGF	hCGa, intact	HVEM	IL-13 R alpha1	IL-17B	IL-17C	IL-17F	IL-17R
8	EpCAM	ErbB2	Erythropoietin R	FSH	HB-EGF	hCGa, intact	HVEM	IL-13 R alpha1	IL-17B	IL-17C	IL-17F	IL-17R
9	Procalcitonin	PSA-free	S-100b	Shh N	Thyroglobulin	Ubiquitin+1	Blank	IC 1	IC 2	IC 3	NEG	NEG
10	Procalcitonin	PSA-free	S-100b	Shh N	Thyroglobulin	Ubiquitin+1	Blank	IC 1	IC 2	IC 3	NEG	NEG

We also offer Custom Human Cytokine Antibody Arrays. You can select the cytokines of interest from the following list and we will produce the customized array at an affordable price. For more information, please visit our website, www.raybiotech.com.

#### **Human Custom Antibody Array List (285 proteins)**

4-18B/TNFRSF9   CNTF   GDNF   IL-18 R alpha   MIP-1 alpha   SCF
Activin A         CRP         GITR Ligand         IL-1ra         MIP-1 delta         SDF-1 alpha           Adiponectin/Acrp30         CTACK/CCL27         GM-CSF         IL-2         MIP-3 alpha         SDF-1 beta           Adipsin/Factor D         CTLA-4         GRO         IL-2 R alpha         MIP-3 beta         sgp130           AFP         CXCL16         GRO-a         IL-2 R beta         MMP-1         Shh N           AgRP(ART)         DAN         Growth Hormom         IL-2 R gamma         MMP-2         Siglec-5           ALCAM         Decorin         HB-EGF         IL-21 R         MMP-3         Siglec-9           Angiogenin         DKK-1         HCC-4/CCL16         IL-22         MMP-7         STNF RII           Angiopoietin-1         DKK-3         hCGa, intact         IL-28A/IFN-lambda         MMP-8         STNT RI           Angiostatin         DPPIV/CD26         HVEM         IL-3         MMP-10         TACE           ANGPTL4         DR6         I-309         IL-31         MMP-13         TECK/CCL26           AR (amphiregulin)         Dtk         ICAM-1         IL-4         MPIF-1         TGF-alpha           AXI         E-Cadherin         ICAM-2         IL-5         MSP a Chain
Adiponectin/Acrp30         CTACK/CCL27         GM-CSF         IL-2         MIP-3 alpha         SDF-1 beta           Adipsin/Factor D         CTLA-4         GRO         IL-2 R alpha         MIP-3 beta         sgp130           AFP         CXCL16         GRO-a         IL-2 R beta         MMP-1         Shh N           AgRP(ART)         DAN         Growth Hormom         IL-2 R gamma         MMP-2         Siglec-5           ALCAM         Decorin         HB-EGF         IL-21 R         MMP-3         Siglec-9           Angiogenin         DKK-1         HCC-4/CCL16         IL-22         MMP-7         sTNF RII           Angiopoietin-1         DKK-3         hCGa, intact         IL-28A/IFN-lambda         MMP-8         sTNT RI           Angiopoietin-2         DKK-4         HGF         IL29/IFN-lambda 1         MMP-9         TACE           Angostatin         DPPIV/CD26         HVEM         IL-31         MMP-10         TARC           ANGPTL4         DR6         I-309         IL-31         MMP-13         TECK/CCL26           AR (amphiregulin)         Dtk         ICAM-1         IL-4         MPIF-1         TGF-alpha           AxI         E-Cadherin         ICAM-2         IL-5         MSP a Chain
Adipsin/Factor D         CTLA-4         GRO         IL-2 R alpha         MIP-3 beta         sgp130           AFP         CXCL16         GRO-a         IL-2 R beta         MMP-1         Shh N           AgRP(ART)         DAN         Growth Hormom         IL-2 R gamma         MMP-2         Siglec-5           ALCAM         Decorin         HB-EGF         IL-21 R         MMP-3         Siglec-9           Angiogenin         DKK-1         HCC-4/CCL16         IL-22         MMP-7         sTNF RII           Angiopoietin-1         DKK-3         hCGa, intact         IL-28A/IFN-lambda         MMP-8         sTNT RI           Angiopoietin-2         DKK-4         HGF         IL29/IFN-lambda 1         MMP-9         TACE           Angiostatin         DPPIV/CD26         HVEM         IL-3         MMP-10         TARC           ANGPTL4         DR6         I-309         IL-31         MMP-13         TECK/CCL26           AR (amphiregulin)         Dtk         ICAM-1         IL-4         MPIF-1         TGF-alpha           AxI         E-Cadherin         ICAM-2         IL-5         MSP a Chain         TGF-beta 1           B7-1(CD80)         EDA-A2         ICAM-3         IL-5 R alpha         NAP-2         TGF-b
AFP         CXCL16         GRO-a         IL-2 R beta         MMP-1         Shh N           AgRP(ART)         DAN         Growth Hormom         IL-2 R gamma         MMP-2         Siglec-5           ALCAM         Decorin         HB-EGF         IL-21 R         MMP-3         Siglec-9           Angiogenin         DKK-1         HCC-4/CCL16         IL-22         MMP-7         sTNF RII           Angiopoietin-1         DKK-3         hCGa, intact         IL-28A/IFN-lambda         MMP-8         sTNT RI           Angiopoietin-2         DKK-4         HGF         IL29/IFN-lambda 1         MMP-9         TACE           Angiostatin         DPPIV/CD26         HVEM         IL-3         MMP-10         TARC           ANGPTL4         DR6         I-309         IL-31         MMP-13         TECK/CCL25           AR (amphiregulin)         Dtk         ICAM-1         IL-4         MPIF-1         TGF-alpha           AxI         E-Cadherin         ICAM-2         IL-5         MSP a Chain         TGF-beta 1           B7-1(CD80)         EDA-A2         ICAM-3         IL-5 R alpha         NAP-2         TGF-beta 2
AgRP(ART)         DAN         Growth Hormom         IL-2 R gamma         MMP-2         Siglec-5           ALCAM         Decorin         HB-EGF         IL-21 R         MMP-3         Siglec-9           Angiogenin         DKK-1         HCC-4/CCL16         IL-22         MMP-7         STNF RII           Angiopoietin-1         DKK-3         hCGa, intact         IL-28A/IFN-lambda         MMP-8         STNT RI           Angiopoietin-2         DKK-4         HGF         IL29/IFN-lambda 1         MMP-9         TACE           Angiostatin         DPPIV/CD26         HVEM         IL-3         MMP-10         TARC           ANGPTL4         DR6         I-309         IL-31         MMP-13         TECK/CCL26           AR (amphiregulin)         Dtk         ICAM-1         IL-4         MPIF-1         TGF-alpha           AxI         E-Cadherin         ICAM-2         IL-5         MSP a Chain         TGF-beta 1           B7-1(CD80)         EDA-A2         ICAM-3         IL-5 R alpha         NAP-2         TGF-beta 2
ALCAM Decorin HB-EGF IL-21 R MMP-3 Siglec-9  Angiogenin DKK-1 HCC-4/CCL16 IL-22 MMP-7 STNF RII  Angiopoietin-1 DKK-3 hCGa, intact IL-28A/IFN-lambda MMP-8 STNT RI  Angiopoietin-2 DKK-4 HGF IL29/IFN-lambda 1 MMP-9 TACE  Angiostatin DPPIV/CD26 HVEM IL-3 MMP-10 TARC  ANGPTL4 DR6 I-309 IL-31 MMP-13 TECK/CCL26  AR (amphiregulin) Dtk ICAM-1 IL-4 MPIF-1 TGF-alpha  AxI E-Cadherin ICAM-2 IL-5 MSP a Chain TGF-beta 1  B7-1(CD80) EDA-A2 ICAM-3 IL-5 R alpha NAP-2 TGF-beta 2
Angiogenin         DKK-1         HCC-4/CCL16         IL-22         MMP-7         sTNF RII           Angiopoietin-1         DKK-3         hCGa, intact         IL-28A/IFN-lambda         MMP-8         sTNT RI           Angiopoietin-2         DKK-4         HGF         IL29/IFN-lambda 1         MMP-9         TACE           Angiostatin         DPPIV/CD26         HVEM         IL-3         MMP-10         TARC           ANGPTL4         DR6         I-309         IL-31         MMP-13         TECK/CCL25           AR (amphiregulin)         Dtk         ICAM-1         IL-4         MPIF-1         TGF-alpha           AxI         E-Cadherin         ICAM-2         IL-5         MSP a Chain         TGF-beta 1           B7-1(CD80)         EDA-A2         ICAM-3         IL-5 R alpha         NAP-2         TGF-beta 2
Angiopoietin-1         DKK-3         hCGa, intact         IL-28A/IFN-lambda         MMP-8         sTNT RI           Angiopoietin-2         DKK-4         HGF         IL29/IFN-lambda 1         MMP-9         TACE           Angiostatin         DPPIV/CD26         HVEM         IL-3         MMP-10         TARC           ANGPTL4         DR6         I-309         IL-31         MMP-13         TECK/CCL25           AR (amphiregulin)         Dtk         ICAM-1         IL-4         MPIF-1         TGF-alpha           AxI         E-Cadherin         ICAM-2         IL-5         MSP a Chain         TGF-beta 1           B7-1(CD80)         EDA-A2         ICAM-3         IL-5 R alpha         NAP-2         TGF-beta 2
Angiopoietin-2         DKK-4         HGF         IL29/IFN-lambda 1         MMP-9         TACE           Angiostatin         DPPIV/CD26         HVEM         IL-3         MMP-10         TARC           ANGPTL4         DR6         I-309         IL-31         MMP-13         TECK/CCL25           AR (amphiregulin)         Dtk         ICAM-1         IL-4         MPIF-1         TGF-alpha           AxI         E-Cadherin         ICAM-2         IL-5         MSP a Chain         TGF-beta 1           B7-1(CD80)         EDA-A2         ICAM-3         IL-5 R alpha         NAP-2         TGF-beta 2
Angiostatin         DPPIV/CD26         HVEM         IL-3         MMP-10         TARC           ANGPTL4         DR6         I-309         IL-31         MMP-13         TECK/CCL25           AR (amphiregulin)         Dtk         ICAM-1         IL-4         MPIF-1         TGF-alpha           AxI         E-Cadherin         ICAM-2         IL-5         MSP a Chain         TGF-beta 1           B7-1(CD80)         EDA-A2         ICAM-3         IL-5 R alpha         NAP-2         TGF-beta 2
ANGPTL4 DR6 I-309 IL-31 MMP-13 TECK/CCL25 AR (amphiregulin) Dtk ICAM-1 IL-4 MPIF-1 TGF-alpha AxI E-Cadherin ICAM-2 IL-5 MSP a Chain TGF-beta 1 B7-1(CD80) EDA-A2 ICAM-3 IL-5 R alpha NAP-2 TGF-beta 2
AR (amphiregulin)         Dtk         ICAM-1         IL-4         MPIF-1         TGF-alpha           AxI         E-Cadherin         ICAM-2         IL-5         MSP a Chain         TGF-beta 1           B7-1(CD80)         EDA-A2         ICAM-3         IL-5 R alpha         NAP-2         TGF-beta 2
AxI         E-Cadherin         ICAM-2         IL-5         MSP a Chain         TGF-beta 1           B7-1(CD80)         EDA-A2         ICAM-3         IL-5 R alpha         NAP-2         TGF-beta 2
B7-1(CD80) EDA-A2 ICAM-3 IL-5 R alpha NAP-2 TGF-beta 2
Rate 2 M FGF   IFN gamma   II .
Indicates in in-Admining lines into information in incompanies in
BCAM EGF R IGFBP-1 IL-6 sR NGF R Thyroglobuli
BCMA/TNFRSF17 EG-VEGF/PK1 IGFBP-2 IL-7 Nidogen-1/Entactin Tie-1
BDNF ENA-78 IGFBP-3 IL-8 NrCAM Tie-2
beta IG-H3 Endoglin IGFBP-4 IL-9 NRG1-beta 1/HRG1-beta 1 TIM-1
Betacellulin (BTC) Endostatin IGFBP-5 IL-9 R NT-3 TIMP-1
bFGF Eotaxin IGFBP-6 Insulin NT-4 TIMP-2
BLC Eotaxin-2 IGF-I IP-10 Oncostatin M TIMP-4
BMP-4 Eotaxin-3 IGF-I sR I-TAC/CXCL11 Osteopontin TNF-alpha
BMP-5 EpCAM/TROP1 IGF-II LAP(TGF-b1) Osteoprotegerin TNF-beta
BMP-6 ErbB2 IL-1 alpha Leptin R PAI-I TPO
BMP-7 ErbB3 IL-1 beta LEPTIN(OB) PARC TRAIL R1
b-NGF Erythropoietin R (EPO R) IL-1 R4/ST2 LH P-Cadherin TRAIL R2
BTC E-Selectin IL-1 sRI LIF PDGF R alpha TRAIL R3
CA125 Fas Ligand IL-1 sRII LIGHT PDGF R beta TRAIL R4
CA15-3 Fas/TNFRSF6 IL-10 LIMPII/SR-B2 PDGF-AA Trappin-2/El.
CA19-9 Fcr RIIB/C IL-10 R alpha Lipocalin-2/NGAL PDGF-AB TREM-1
Carbonic Anhydrase IX(CA9) Ferritin IL-10 R beta L-Selectin PDGF-BB TROY
Cardiotrophin-1 (CT-1) FGF-4 IL-11 Lymphotactin PECAM-1 TSH
Cathepsin S FGF-6 IL-12 p40 LYVE-1 Platelet Factor 4 TSLP
CCL14a/HCC-1 FGF-7 IL-12 p70 Marapsin/Pancreasin PIGF u PAR
CCL21/6ckine FGF-9 IL-13 MCP-1 Procalcitonin/Calcitonin Ubiquitin+1
CCL28/VIC FLRG IL-13 Ra1 MCP-2 Prolactin VCAM-1
CD14 Fit-3 Ligand IL-13 Ra2 MCP-3 PSA-free VE-Cadherin
CD23/Fc epsilon RII Follistatin IL-15 MCP-4 PSA-total VEGF
CD27 Fractalkine IL-16 MCSF P-selectin VEGF R2
,
CD30 FSH IL-17 M-CSF R RAGE VEGF R3
CD30         FSH         IL-17         M-CSF R         RAGE         VEGF R3           CD40         Furin         IL-17B         MDC         RANK         VEGF-C
CD40 Furin IL-17B MDC RANK VEGF-C
CD40         Furin         IL-17B         MDC         RANK         VEGF-C           CD40 Ligand         Galectin-7         IL-17C         MICA         RANTES         VEGF-D

RayBiotech, Inc., the protein array pioneer company, strives to research and develop new products to meet demands of the biomedical community. RayBio's patent-pending technology allows detection of over 180 cytokines, chemokines and other proteins in a single experiment. Our format is simple, sensitive, reliable and cost effective. Products include: Cytokine Arrays, Chemokine Arrays, ELISA kits, Phosphotyrosine kits, Recombinant Proteins, Antibodies, and custom services.

#### 1. Antibody arrays

Cytokine antibody array

Human cytokine antibody arrays

Mouse cytokine antibody arrays

Rat cytokine antibody arrays

Pathway- or disease-focused antibody arrays

Inflammation antibody array

Angiogensis antibody array

Chemokine antibody array

Growth factor antibody array

MMP antibody array

Atherosclerosis antibody array

Quantibody arrays for quantitative measurement of cytokine and other protein concentration

Phosphorylation antibody arrays

Biotin label-based antibody arrays for high density antibody arrays.

Antibody analysis tool, software

- 2. ELISA
- 3. Cell-based phosphorylation assay
- 4. Custom antibody arrays
- 5. Antibody
- 6. Recombinant protein
- 7. Peptide
- 8. Protein arrays
- 9. EIA

RayBiotech also provides excellent custom service:

- 1. Antibody arrays
- 2. Protein arrays
- 3. Peptide synthesis

- 4. Production of recombinant protein and antibody
- 5. Peptide arrays
- 6. Phosphorylation arrays
- 7. ELISA
- 8. EIA

Just simply send your samples and we will do the assay for you.

#### Technology transfer program

Have you developed technologies or reagents interested to the scientific and research community? RayBiotech can help you commercialize your technologies, reagents and dream.

# VI. Troubleshooting guide

Problem	Cause	Recommendation
Weak signal	Inadequate detection	Check laser power and PMT parameters
	Inadequate reagent volumes or improper dilution	Check pipetters and ensure correct preparation
	Short incubation times	Ensure sufficient incubation Time and change sample incubation step to overnight
	Too low protein concentration in sample	Don't make too low dilution Or concentrate sample
	Improper storage of kit	Store kit at suggested temperature
High background	Excess of biotinylated antibodies	Make sure correct amount of antibodies
	Excess of streptavidin	Make sure correct amount of streptavidin
	Inadequate detection	Check laser power And PMT parameters
	dust	Work in clean environment
	Insufficient wash	Increase wash time and use more wash buffer
Uneven signal	Bubbles formed during incubation	Avoid bubble formation during incubation
	Arrays are not completed Covered by reagent	Completely cover arrays with solution

#### VII. Selected References Using RayBiotech Products:

Classification and prediction of clinical Alzheimer's diagnosis based on plasma signaling proteins. Ray S, Britschgi M, Herbert C, Takeda-Uchimura Y, et al. **Nature Med**. 2007; 13(11):1359-1362.

Human AB Serum and Thrombin-Activated Platelet-Rich Plasma Are Suitable Alternatives to Fetal Calf Serum for the Expansion of Mesenchymal Stem Cells from Adipose Tissue. Kocaoemer A, Kern S, Kluter H, Bieback K. **Stem Cells**. 2007; 25:1270-1278

ATP Binding by Monarch-1/NLRP12 Is Critical for Its Inhibitory Function. Ye Z, Lich JD, Moore CB, Duncan JA, Williams KL, Ting JP-Y. **Mol Cell Biol**. 2008;28:1841-1850..

Systemic Endocrine Instigation of Indolent Tumor Growth Requires Osteopontin. McAllister SS, Gifford AM, Greiner AL, Kelleher SP, Saelzer MP, et al. **Cell**. 2008;133:994-1005.

Chemokine Signaling via the CXCR2 Receptor Reinforces Senescence. Acosta JC, O'Loghlen A, Banito A, Guijarro MV, Augert A, et al. **Cell**. 2008;133: 1006-1018.

HIV-1-mediated apoptosis of neuronal cells: Proximal molecular mechanisms of HIV-1-induced encephalopathy. Yan Xu, Joseph Kulkoshy, Roger J. Pomerantz. **PNAS**. 2004 May 4, 2004 Vol. 101 No. 18.

Veto-like activity of mesenchymal stem cells: functional discrimination between cellular responses to alloantigens and recall antigens. Rameshwar P. **Journal of Immunology**. 2003 Oct 1;171(7):3426-34.

Monocyte Chemotactic Protein-1 Mediates Prostate Cancer-Induced Bone Resorption. Lu Y, Cai Z, Xiao G, Keller ET, et al. **Cancer Res**. 2007; 67(8):3646-3653.

Proteomic Characterization of the Interstitial Fluid Perfusing the Breast Tumor Microenvironment. A Novel Resource for Biomarker and Therapeutic Target Discovery. Julio E. Celis, Pavel Gromov, Teresa Cabezón, José M. A. Moreira, Noona Ambartsumian, Kerstin Sandelin, Fritz Rank, and Irina Gromova. **Molecular Cellular Proteomics**. April 2004; 11(3):328-39.

Antibody array-generated profiles of cytokine release from THP-1 leukemic monocytes exposed to different amphotericin B formulations. Turtinen LW, Prall DN, Bremer LA, Nauss RE, Hartsel SC. **Antimicrobial Agents Chemotherapy**. 2004 Feb;48(2):396-403.

Reduced T-cell and dendritic cell function is related to cyclooxygenase-2 overexpression and protaglandin e(2) secretion in patients with breast cancer. Pockaj BA, Basu GD. **Annals of Surgical Oncology**. 3:327-344, 2004.

Inhibition of macrophage migration inhibitory factor decreases proliferation and cytokine expression in bladder cancer cells. Katherine L Meyer-Siegler, **BMC Cancer**. 2004, 4:34.

Airway cytokine expression measured by means of protein array in exhaled breath condensate: Correlation with physiologic properties in asthmatic patients. Matsunaga K, Yanagisawa S, Ichikawa T, Ueshima K, et al. **J Allergy Clin Immunol**. 2006; 188:84-90.

In Vivo Proteomic Analysis of Cytokine Expression in Laser Capture-Microdissected Urothelial Cells of Obstructed Ureteropelvic Junction Procured by Laparoscopic Dismembered Pyeloplasty. **Journal of Endourology**. 2003 June; Volume:17 Number:5 Page:333-336.

#### Note:

RayBio<sup>®</sup> is the trademark of RayBiotech, Inc.

Cytokine protein arrays are RayBiotech patent-pending technology.

This product is intended for research only and is not to be used for clinical diagnosis. Our produces may not be resold, modified for resale, or used to manufacture commercial products without written approval by RayBiotech, Inc.

Under no circumstances shall RayBiotech be liable for any damages arising out of the use of the materials.

Products are guaranteed for three months from the date of purchase when handled and stored properly. In the event of any defect in quality or merchantability, RayBiotech's liability to buyer for any claim relating to products shall be limited to replacement or refund of the purchase price.

HiLytePlus™ is a trademark of AnaSpec, Inc.

This product is for research use only.



©2009 RayBiotech, Inc.