## University of Iowa NMR Central Research Facility World Wide Web Scheduling User's Manual

Prepared 27 June, 2001

University of Iowa High Field NMR Facility Iowa City, Iowa 52242 (319) 335-1332

This manual is intended for use by the University of Iowa High Field NMR Facility's users. Unauthorized use, access, or tampering is strictly forbidden. Use of this manual and the internet pages described within constitutes acceptance of all of the Facility's polices. The University and the Facility accepts no responsibility for lost data or damage to any hardware that results from misuse of the internet pages described herein.

The NMR Central Research Facility accepts reservations for time via its homepage on the World Wide Web. In order to reserve time, you must be checked out on the instrument, and have an account. See John Snyder or Dr. Donald Stec if you are checked out but do not have an account. If you are not checked out, see John Snyder for the Facility's polices regarding instrument check out.

Assuming you have an account, begin by starting any web browser and opening the location:

http://nmr.chem.uiowa.edu

Your browser should pull up the Facility's homepage:



Figure 1: The NMR Facility's homepage

Click on the word <u>schedule</u> on the left of the screen. This will bring up the scheduling homepage. Among other things, you may sign on and reserve time from this page:



Figure 2: The scheduling page.

Click: "<u>Click Here to Reserve Time</u>". From here, you will see the login page:



Figure 3: Login page

Here, enter your login name and password in the spaces provided. Then click Login. This will bring up the instrument and day selection screen.



Figure 4: Instrument and day selection screen.

If you do not get this screen, note the error, and call John Snyder. From the pull-down menus, select the instrument and day that you need, then click Step Two.

逝い	Iniver	sity a	fIow	a NMI	t Central	Resear	ch Facility	- Netscape							_ 🗆 ×
Ele	Edit	View	<u>G</u> o	⊆orn	nunicator	Help									
	aci Baci	. F	ر orwar	rd F	3. ieload	A Home	arch 🌽	Metscape	it Brint	Security SH	). hop	Stop			N
	¥.	Bookn	arks	4	Netsite:	http://nmr	chem.uiow	a.edu/					· 🗇	What	's Related
University of Iowa												litv			
Home											_				
	Schedule Timesheat for the DPV300 on												-		
-	Timesneet for the DPA500 on														
	Staff Saturday														
	P	olic	y												
	nsti	um	ent	ts					1		-1				
								Time	Status	Action					
	IVIa	inu	ais					800	open	C reserve	-1				
	F	orm	s					830	open	C reserve	-1-				
F		ink	•					900	open	C reserve					
		IIIK	•					930	open	C reserve					
								1000	walk	Not Availat	ble				
								1010	walk	Not Availat	ble				
								1020	walk	Not Availat	ble				
								1030	walk	Not Availa	ble				
								1040	walk	Not Availat	ble				
								1050	walk	Not Availat	ole				
								1100	walk	Not Availat	ble				•
							[Herae]	[Schedule] [Stat	f] [Policy] []	nstruments] [Mun	utals] [Fen	ns) (Lir	de ]		
8	-0				Documen	t: Done					- <b>**</b>	9 <u>85</u>	d۵	<b></b>	1

Figure 5: Time reservation screen.

If open appears in the Action column, the instrument is available, and you may reserve time, by clicking on the "radio button" next to the word reserve in the Action column. When all the time you need is reserved, click the Execute button at the bottom of the sheet:



Figure 6: New timesheet.

That's it. You're done. The time you entered on the previous page is now listed in the new timesheet. From here, you can quit your browser, or continue to browse other pages. If you need to cancel time, log on as usual, and any time that you have reserved will have the word "unreserve" next to it in the Action column. Click its button as if you were reserving time, and click the Execute button. The time will be canceled.

💥 University of Iowa NMR Central Research Facility - Netscape															
Ele	Edit	View	Go	⊆ommu	nicator	Help									
-	Back		ionward	i Rel	) pad	And Home	🥭 Search	Netscape	i d 🕉 Print	ing Security	🔕. Shop	Stop		Ν	
1	<b>(</b>	Bookm	arks	🧶 N	Netsite: http://mm.chem.uiowa.edu/								🗊 " What'	s Related	
University of Iowa NMR Central Research Facilit													lty		
	Home Timesheat for the DPV300 on												-		
													-		
	Timesneet for the DFA500 on														
	Staff Saturday														
	Р	olic	у	_											
	nstr	um	ent	s											
H	Ma	nur	ale					Time	Status	Actio	n				
	IVIC	nua	115					800	luci	• unrese	rve				
	F	orm	S					830	luci	• unrese	rve				
	1	ink						900	open	C reserve					
			<u> </u>					930	open	C reserve					
L								1000	walk	Not Avail	able				
L								1010	walk	Not Avail	able				
L								1020	walk	Not Avail	able				
L								1030	walk	Not Avail	able				
								1040	walk	Not Avail	able				
								1050	walk	Not Avail	able				
								1100	walk	Not Avail	able			-	
[Henne] [Schednike] [Staff] [Policy] [Instruments] [Manuals] [Ferrar] [Limbs]															
2	<b>0</b> -1			D	ocumen	t: Done					- 💥	, 🤒 🥶	P 👔	1.	

If you encounter any difficulties with this service, please feel free to contact the Facility at (319) 335-1332.