



Genomic DNA from blood

User manual

NucleoMag[®] Blood 3 mL

March 2012/Rev.02

Table of contents

1	Components	4
1.1	Kit contents	4
1.2	Material to be supplied by the user	4
2	Product description	5
2.1	The basic principle	5
2.2	Kit specifications	5
2.3	Elution procedures	6
3	Storage conditions and preparation of working solutions	7
4	Safety instructions	8
4.1	Risk and safety phrases	8
4.2	GHS classification	9
5	Protocol for the isolation of genomic DNA from 3 mL blood samples using KingFisher® Flex 24	12
6	Appendix	14
6.1	Troubleshooting	14
6.3	Product use restriction/warranty	16

1 Components

1.1 Kit contents

NucleoMag® Blood 3 mL	
REF	1x 96 preps 744502.1
NucleoMag® B-Beads	18 mL
Lysis Buffer MBL1	125 mL
Binding Buffer MBL2	550 mL
Wash Buffer MBL3	1 000 mL
Wash Buffer MBL4	500 mL
Elution Buffer MBL5*	125 mL
Proteinase K, lyophilized**	12 x 75 mg
Proteinase Buffer PB	2 x 35 mL
User manual	1

1.2 Material to be supplied by the user

- Magnetic separator
KingFisher® Flex 24 instrument
- Separation plates, elution plates
KingFisher® 24 deep well plates
- KingFisher® 24 well tip comb
- 80 % ethanol (for the washing step)

* Elution Buffer MBL5: 5 mM Tris, pH 8.5

**For preparation of working solutions and storage conditions see section 3.

2 Product description

2.1 The basic principle

The **NucleoMag® Blood 3 mL** procedure is based on reversible adsorption of nucleic acids to paramagnetic beads under appropriate buffer conditions. Whole blood is lysed with Lysis Buffer MBL1 and Proteinase K. Following lysis incubation, magnetic beads are added and binding conditions under which the DNA binds to the magnetic beads are adjusted by addition of Binding Buffer MBL2. After magnetic separation and removal of supernatant, the paramagnetic beads are washed three times to remove contaminants and salt. There is no need for a drying step as ethanol from previous wash steps is removed by Wash Buffer MBL4. Finally, highly purified DNA is eluted with low-salt Elution Buffer MBL5 and can directly be used for downstream applications. **NucleoMag® Blood 3 mL** is recommended for use on KingFisher® Flex 24 instrument.

2.2 Kit specifications

The **NucleoMag® Blood 3 mL** kit is made for isolation of genomic DNA from blood samples. This kit provides reagents and magnetic beads for isolation of genomic DNA from 96 samples of up to 3 mL. The purified DNA can be used directly as template for PCR, blotting, or any kind of enzymatic reactions.

The kit provides reagents for the purification of up to 100–130 µg of pure genomic DNA from 3 mL whole blood with an A_{260}/A_{280} ratio ≥ 1.6 –1.9.

Fresh, frozen, or blood treated either with EDTA or citrate can be used.

NucleoMag® Blood 3 mL kit can be processed completely at room temperature. Elution at 55 °C will increase the yield by about 15–20 %.

NucleoMag® B-Beads are highly reactive, superparamagnetic beads with a high binding capacity.

NucleoMag® Blood 3 mL kit has been developed for use with ThermoFisher's KingFisher® Flex 24 instrument. A script is available on request from MACHEREY-NAGEL. The maximum sample volume of 3 mL is splitted into two aliquots of 1.5 mL each.

For processing smaller blood sample volumes, use of liquid handling robots other than the KingFisher® Flex 24 or manual extraction, please inquire with MN technical support for details.

For smaller blood sample volumes, MN offers the **NucleoMag® Blood 200 µL** kit (see ordering information).

2.3 Elution procedures

Purified genomic DNA can be eluted directly with the supplied Elution Buffer MBL5. Elution can be carried out in a volume of > 1 mL. Smaller elution buffer volumes may result in incomplete bead separation. For efficient elution, the magnetic bead pellet should be resuspended completely in the elution buffer.

3 Storage conditions and preparation of working solutions

Attention: Buffers MBL1, MBL2, and MBL3 contain chaotropic salt! Wear gloves and goggles!

CAUTION: Buffer MBL1 contains guanidine hydrochloride which can form highly reactive compounds when combined with bleach (sodium hypochlorite). DO NOT add bleach or acidic solutions directly to the sample-preparation waste.

- All components of the **NucleoMag® Blood 3 mL** kit should be stored at room temperature (18–25 °C) and are stable for up to one year.
- All buffers are delivered ready-to-use.

Before starting **NucleoMag® Blood 3 mL** protocol prepare the following:

- Add the indicated volume of Proteinase Buffer PB to dissolve lyophilized Proteinase K (see table below). Proteinase K solution is stable at -20 °C for up to 6 months.

NucleoMag® Blood 3 mL	
REF	1 x 96 preps 744502.1
Proteinase K	Add 3.75 mL Proteinase Buffer PB to each vial

4 Safety instructions

The following components of the **NucleoMag® Blood 3 mL** kits contain hazardous contents.

Wear gloves and goggles and follow the safety instructions given in this section.

4.1 Risk and safety phrases

Component	Hazard contents	Hazard symbol	Risk phrases	Safety phrases
<i>Inhalt</i>	<i>Gefahrstoff</i>	<i>Gefahrstoff-symbol</i>	<i>R-Sätze</i>	<i>S-Sätze</i>
MBL1	Guanidine hydrochloride 50–66 % <i>Guanidinhydrochlorid 50–66 %</i>	✘ Xi*	R 22-36/38	S 26-37/39
MBL2	Sodium perchlorate 20–40 % + ethanol 35–55 % <i>Natriumperchlorat 20–40 % + Ethanol 35–55 %</i>	*	R 10-22	S 13-16
MBL3	Sodium perchlorate 5–20 % + ethanol 20–35 % <i>Natriumperchlorat 5–20 % + Ethanol 20–35 %</i>	*	R 10	S 16
Proteinase K	Proteinase K, lyophilized <i>Proteinase K, lyophilisiert</i>	✘ Xn	R 36/37/38-42	S 22-24-26-36/37

Risk phrases

R 10	Flammable. <i>Entzündlich.</i>
R 22	Harmful if swallowed. <i>Gesundheitsschädlich bei Verschlucken.</i>
R 36/37/38	Irritating to eyes, respiratory system and skin. <i>Reizt die Augen, Atmungsorgane und die Haut.</i>
R 36/38	Irritating to eyes and skin. <i>Reizt die Augen und die Haut.</i>
R 42	May cause sensitization by inhalation. <i>Sensibilisierung durch Einatmen und Hautkontaktmöglich.</i>

* Hazard labeling not necessary if quantity per bottle below 125 g or mL (certificate of exemption according to 67/548/EEC Art. 25, 1999/45/EC Art. 12 and German GefStoffV § 20 (3) and TRGS 200 7.1). For further information see Material Safety Data Sheet.





Safety phrases

- S 13 Keep away from food, drink and animal foodstuffs.
Von Nahrungsmitteln, Getränken und Futtermitteln fernhalten.
- S 16 Keep away from sources of ignition – No smoking!
Von Zündquellen fernhalten – Nicht rauchen!
- S 22 Do not breathe dust.
Staub nicht einatmen.
- S 24 Avoid contact with the skin.
Berührung mit der Haut vermeiden.
- S 26 In case of contact with eyes, rinse immediately with plenty of water and seek medical advice.
Bei Berührung mit den Augen sofort gründlich mit Wasser abspülen und Arzt konsultieren.
- S 36/37 Wear suitable protective clothing and gloves.
Bei der Arbeit geeignete Schutzhandschuhe und Schutzbrille / Gesichtsschutz tragen.
- S 37/39 Wear suitable gloves and eye / face protection.
Bei der Arbeit geeignete Schutzhandschuhe und Schutzkleidung tragen.

4.2 GHS classification

Only harmful features do not need to be labeled with H and P phrases until 125 mL or 125 g.

Mindergefährliche Eigenschaften müssen bis 125 mL oder 125 g nicht mit H- und P-Sätzen gekennzeichnet werden.

Component	Hazard contents	GHS symbol	Hazard phrases	Precaution phrases
<i>Inhalt</i>	<i>Gefahrstoff</i>	<i>GHS Symbol</i>	<i>H-Sätze</i>	<i>P-Sätze</i>
MBL1	Guanidine hydrochloride 50–66 % <i>Guanidinhydrochlorid 50–66 %</i>	 Danger <i>Gefahr</i>	302, 315, 319	280, 301+312, 302+352, 305+351+338, 330, 332+313, 337+313
MBL2	Sodium perchlorate 20–40 % + ethanol 35–55 % <i>Natriumperchlorat 20–40 % + Ethanol 35–55 %</i>	 Warning <i>Achtung</i>	226, 302	210, 233, 301+312, 330, 403+235
MBL3	Sodium perchlorate 5–20 % + ethanol 20–35 % <i>Natriumperchlorat 5–20 % + Ethanol 20–35 %</i>	 Warning <i>Achtung</i>	226	210, 233, 403+235
Proteinase K	Proteinase K, lyophilized <i>Proteinase K, lyophilisiert</i>	 Danger <i>Gefahr</i>	315, 319, 334, 335	261, 280, 302+352, 304+340, 305+351+338, 312, 332+313, 337+313, 342+311, 403+233

Hazard phrases

- H 226 Flammable liquid and vapour.
Flüssigkeit und Dampf entzündbar.
- H 302 Harmful if swallowed.
Gesundheitsschädlich bei Verschlucken.
- H 315 Causes skin irritation.
Verursacht Hautreizungen.
- H 319 Causes serious eye irritation.
Verursacht schwere Augenreizung.
- H 334 May cause allergy or asthma symptoms or breathing difficulties if inhaled.
Kann bei Einatmen Allergie, asthmaartige Symptome oder Atembeschwerden verursachen.
- H 335 May cause respiratory irritation.
Kann die Atemwege reizen.

Precaution phrases

- P 210 Keep away from heat/sparks/open flames/hot surfaces – No smoking.
Von Hitze/Funken/offener Flamme/heißen Oberflächen fernhalten. Nicht rauchen.
- P 233 Keep container tightly closed.
Behälter dicht verschlossen halten.
- P 261 Avoid breathing dust.
Einatmen von Staub vermeiden.
- P 280 Wear protective gloves / eye protection.
Schutzhandschuhe/Augenschutz tragen.
- P 301+312 IF SWALLOWED: Call a POISON CENTER or doctor / physician if you feel unwell.
BEI VERSCHLUCKEN: Bei Unwohlsein GIFTINFORMATIONSZENTRUM oder Arzt anrufen.
- P 302+352 IF ON SKIN: Wash with plenty of soap and water.
BEI KONTAKT MIT DER HAUT: Mit viel Wasser und Seife waschen.
- P 304+340 IF INHALED: Remove victim to fresh air and keep at rest in a position comfortable for breathing.
BEI EINATMEN: An die frische Luft bringen und in einer Position ruhigstellen, die das Atmen erleichtert.
- P 305+351+338 IF IN EYES: Rinse continuously with water for several minutes. Remove contact lenses if present and easy to do – continue rinsing
BEI KONTAKT MIT DEN AUGEN: Einige Minuten lang behutsam mit Wasser spülen. Vorhandene Kontaktlinsen nach Möglichkeit entfernen. Weiter spülen.
- P 312 Call a POISON CENTER or doctor / physician if you feel unwell.
Bei Unwohlsein GIFTINFORMATIONSZENTRUM oder Arzt anrufen.
- P 330 Rinse mouth.
Mund ausspülen.
- P 332+313 IF skin irritation occurs: Get medical advice / attention.
Bei Hautreizung: Ärztlichen Rat einholen / ärztliche Hilfe hinzuziehen.
- P 337+313 Get medical advice / attention.
Bei anhaltender Augenreizung: Ärztliche Rat einholen / ärztliche Hilfe hinzuziehen.
- P 342+311 If experiencing respiratory symptoms: Call a POISON CENTER or doctor / physician.
Bei Symptomen der Atemwege: Giftinformationszentrum oder Arzt anrufen.
- P 403+233 Store in a well ventilated place. Keep cool.
Behälter dicht verschlossen an einem gut belüfteten Ort aufbewahren.

Precaution phrases

P 403+235 Store in a well ventilated place. Keep cool.
Kühl an einem gut belüfteten Ort aufbewahren.

For further information please see Material Safety Data Sheets (www.mn-net.com).

Weiterführende Informationen finden Sie in den Sicherheitsdatenblättern (www.mn-net.com).

5 Protocol for the isolation of genomic DNA from 3 mL blood samples using KingFisher® Flex 24

The script necessary to run the **NucleoMag® Blood 3 mL** kit on the KingFisher® Flex 24 is available through MN technical support.

Before starting the preparation:

- Check if Proteinase K was prepared according to section 3.
-

1 Lyse sample

Prepare KingFisher® 24 deep well plate with buffers (label deep-well blocks before use).

Wash and elution buffers

Fill **1 mL Elution Buffer MBL5** to each well of an empty Thermo KingFisher® 24 deep well plate.

Fill **4.8 mL Wash Buffer MBL4** to each well of an empty Thermo KingFisher® 24 deep well plate.

Fill **4.8 mL 80 % ethanol** to each well of an empty Thermo KingFisher® 24 deep well plate.

Fill **4.8 mL Wash Buffer MBL3** to each well of an empty Thermo KingFisher® 24 deep well plate.

Fill **4.8 mL Wash Buffer MBL3** to each well of a second empty Thermo KingFisher® 24 deep well plate.

Fill **150 µL of Proteinase K working solution** to each well of the two lysis plates (Thermo KingFisher® 24 deep-well plates).

Samples

Please note that 3 mL blood samples have to be split and distributed into two plates (1.5 mL for each plate)!

Fill **1.5 mL blood sample** to a well of the lysis plate (Thermo KingFisher® 24 deep well plate with 150 µL Proteinase K per well). Fill **1.5 mL blood** of the same sample to the well at the same position of the second lysis plate.

Make sure that one sample is distributed into the same position of each deep well plate (e.g., sample 1 to position A1 of lysis plate 1 and position A1 of lysis plate 2; sample 2 to position A2 of lysis plate 1 and position A2 of lysis plate 2, etc.)

After adding the samples add **575 µL Buffer MBL1** to each well of the two lysis plates.

2 Start isolation on King Fisher® Flex 24 instrument

Start method “NucleoMag®_Blood_3mL” (method is available from MN on request).

Insert plates as indicated on the KingFisher® instrument display.

Method starts with a mixing step (sample lysis) after setting up the last plate to the instrument.

After mixing steps for lysis (approx. 10 min) the instrument will ask for addition of Buffer MBL2 and NucleoMag® B-Beads.

3 Addition of Binding Buffer MBL2 and NucleoMag® B-Beads to lysis plate 1

Add **2.3 mL Buffer MBL2** and **150 µL NucleoMag® B-Beads** to each well of the lysis plate 1.

Mix up NucleoMag® B-Beads before use.

Return lysis plate 1 to the instrument and continue.

4 Addition of Binding Buffer MBL2 to lysis plate 2

Add **2.3 mL Buffer MBL2** to each well of the lysis plate 2.

Return lysis plate 2 to the instrument and continue.

All further steps are now processed without further user interaction.

5 Remove eluted DNA

The instrument stops after the final elution step. Follow the instructions on instrument display and unload the plates from the instrument.

Purified DNA should be centrifuged before UV measurement!

6 Appendix

6.1 Troubleshooting

Problem	Possible cause and suggestions
Poor DNA yield	<i>Elution buffer volume insufficient</i> <ul style="list-style-type: none">• Beads pellet must be covered completely with elution buffer.
	<i>Insufficient performance of elution buffer during elution step</i> <ul style="list-style-type: none">• Remove residual buffers during the separation steps completely. Remaining buffers decrease efficiency of subsequent washing and elution steps.
	<i>Beads dried out</i> <ul style="list-style-type: none">• Do not let the beads dry as this might result in lower elution efficiency.
	<i>Partial elution in Wash Buffer MBL4 already</i> <ul style="list-style-type: none">• Do not resuspend beads in Buffer MBL4 and do not incubate beads in this buffer for more than 2 min, as this buffer is water-based and might elute the DNA already.
	<i>Incubation after dispensing beads to lysate</i> <ul style="list-style-type: none">• Mix immediately after dispensing NucleoMag® B-Beads and Binding Buffer MBL2 to the lysate.
Low purity	<i>Poor blood quality</i> <ul style="list-style-type: none">• Be sure that no blood clots are transferred to the lysis plates. Blood can be stored at 2–8 °C for two weeks. Freeze samples if stored for longer periods.
	<i>Incomplete magnetic bead separation</i> <ul style="list-style-type: none">• High amounts of eluted DNA increase the viscosity and prevent the beads from being attracted completely to the magnets. Increase elution buffer volume.

Problem	Possible cause and suggestions
Suboptimal performance of DNA in downstream applications	<p><i>Carry-over of ethanol from ethanol wash step</i></p> <ul style="list-style-type: none"> • Be sure to remove all of the ethanol from the ethanol wash step. Carry-over of ethanol may interfere with downstream applications. Typically washing the beads in Buffer MBL4 is sufficient to remove ethanol. However, if necessary include a 10 min airdrying step following the Buffer MBL4 wash step.
	<p><i>Low purity</i></p> <ul style="list-style-type: none"> • See above
Carry-over of beads	<p><i>Time for magnetic separation too short</i></p> <ul style="list-style-type: none"> • Increase separation time to allow the beads to be completely attracted to the magnets.
	<p><i>Incomplete magnetic bead separation</i></p> <ul style="list-style-type: none"> • High amounts of eluted DNA increase the viscosity and prevent the beads from being attracted completely to the magnets. Increase elution buffer volume.
Cross contamination	<p><i>Overfilling of wells from the 24-well separation plate</i></p> <ul style="list-style-type: none"> • Do not overfill the wells of the separation plates to avoid cross contamination by splashing.

6.2 Ordering information

Product	REF	Pack of
NucleoMag® Blood 3 mL	744502.1	1 x 96 preps
NucleoMag® Blood 200 µL	744501.1	1 x 96 preps
	744501.4	4 x 96 preps

Visit www.mn-net.com for more detailed product information.

6.3 Product use restriction/warranty

NucleoMag® Blood 3 mL kit components are intended, developed, designed, and sold FOR RESEARCH PURPOSES ONLY, except, however, any other function of the product being expressly described in original MACHEREY-NAGEL product leaflets.

MACHEREY-NAGEL products are intended for GENERAL LABORATORY USE ONLY! MACHEREY-NAGEL products are suited for QUALIFIED PERSONNEL ONLY! MACHEREY-NAGEL products shall in any event only be used wearing adequate PROTECTIVE CLOTHING. For detailed information please refer to the respective Material Safety Data Sheet of the product! MACHEREY-NAGEL products shall exclusively be used in an ADEQUATE TEST ENVIRONMENT. MACHEREY-NAGEL does not assume any responsibility for damages due to improper application of our products in other fields of application. Application on the human body is STRICTLY FORBIDDEN. The respective user is liable for any and all damages resulting from such application.

DNA/RNA/PROTEIN purification products of MACHEREY-NAGEL are suitable for *IN VITRO*-USES ONLY!

ONLY MACHEREY-NAGEL products specially labeled as IVD are also suitable for *IN VITRO*-diagnostic use. Please pay attention to the package of the product. *IN VITRO*-diagnostic products are expressly marked as IVD on the packaging.

IF THERE IS NO IVD SIGN, THE PRODUCT SHALL NOT BE SUITABLE FOR *IN VITRO*-DIAGNOSTIC USE!

ALL OTHER PRODUCTS NOT LABELED AS IVD ARE NOT SUITED FOR ANY CLINICAL USE (INCLUDING, BUT NOT LIMITED TO DIAGNOSTIC, THERAPEUTIC AND/OR PROGNOSTIC USE).

No claim or representations is intended for its use to identify any specific organism or for clinical use (included, but not limited to diagnostic, prognostic, therapeutic, or blood banking). It is rather in the responsibility of the user or - in any case of resale of the products - in the responsibility of the reseller to inspect and assure the use of the DNA/RNA/protein purification products of MACHEREY-NAGEL for a well-defined and specific application.

MACHEREY-NAGEL shall only be responsible for the product specifications and the performance range of MN products according to the specifications of in-house quality control, product documentation and marketing material.

This MACHEREY-NAGEL product is shipped with documentation stating specifications and other technical information. MACHEREY-NAGEL warrants to meet the stated specifications. MACHEREY-NAGEL's sole obligation and the customer's sole remedy is limited to replacement of products free of charge in the event products fail to perform as warranted. Supplementary reference is made to the general business terms and conditions of MACHEREY-NAGEL, which are printed on the price list. Please contact us if you wish to get an extra copy.

There is no warranty for and MACHEREY-NAGEL is not liable for damages or defects arising in shipping and handling (transport insurance for customers excluded), or out of accident or improper or abnormal use of this product; defects in products or

components not manufactured by MACHEREY-NAGEL, or damages resulting from such non-MACHEREY-NAGEL components or products.

MACHEREY-NAGEL makes no other warranty of any kind whatsoever, and SPECIFICALLY DISCLAIMS AND EXCLUDES ALL OTHER WARRANTIES OF ANY KIND OR NATURE WHATSOEVER, DIRECTLY OR INDIRECTLY, EXPRESS OR IMPLIED, INCLUDING, WITHOUT LIMITATION, AS TO THE SUITABILITY, REPRODUCTIVITY, DURABILITY, FITNESS FOR A PARTICULAR PURPOSE OR USE, MERCHANTABILITY, CONDITION, OR ANY OTHER MATTER WITH RESPECT TO MACHEREY-NAGEL PRODUCTS.

In no event shall MACHEREY-NAGEL be liable for claims for any other damages, whether direct, indirect, incidental, compensatory, foreseeable, consequential, or special (including but not limited to loss of use, revenue or profit), whether based upon warranty, contract, tort (including negligence) or strict liability arising in connection with the sale or the failure of MACHEREY-NAGEL products to perform in accordance with the stated specifications. This warranty is exclusive and MACHEREY-NAGEL makes no other warranty expressed or implied.

The warranty provided herein and the data, specifications and descriptions of this MACHEREY-NAGEL product appearing in MACHEREY-NAGEL published catalogues and product literature are MACHEREY-NAGEL's sole representations concerning the product and warranty. No other statements or representations, written or oral, by MACHEREY-NAGEL's employees, agent or representatives, except written statements signed by a duly authorized officer of MACHEREY-NAGEL are authorized; they should not be relied upon by the customer and are not a part of the contract of sale or of this warranty.

Product claims are subject to change. Therefore please contact our Technical Service Team for the most up-to-date information on MACHEREY-NAGEL products. You may also contact your local distributor for general scientific information. Applications mentioned in MACHEREY-NAGEL literature are provided for informational purposes only. MACHEREY-NAGEL does not warrant that all applications have been tested in MACHEREY-NAGEL laboratories using MACHEREY-NAGEL products. MACHEREY-NAGEL does not warrant the correctness of any of those applications.

Last updated: 07/2010, Rev. 03

Please contact:
MACHEREY-NAGEL GmbH & Co. KG
Tel.: +49 24 21 969-270
tech-bio@mn-net.com

Trademarks:

KingFisher® is a registered trademark of Thermo Fisher Scientific

NucleoMag® is a registered trademark of MACHEREY-NAGEL GmbH & Co. KG

All used names and denotations can be brands, trademarks, or registered labels of their respective owner – also if they are not special denotation. To mention products and brands is only a kind of information (i.e., it does not offend against trademarks and brands and can not be seen as a kind of recommendation or assessment). Regarding these products or services we can not grant any guarantees regarding selection, efficiency, or operation.