

**Air Jet Sieve
SV016**

Impact Test Equipment Ltd
www.impact-test.co.uk & www.impact-test.com

User Guide
User Guide
User Guide

**Impact Test Equipment Ltd.
Building 21 Stevenston Ind. Est.
Stevenston
Ayrshire
KA20 3LR**

T: 01294 602626

F: 01294 461168

E: sales@impact-test.co.uk

Test Equipment Web Site

www.impact-test.co.uk

Test Sieves & Accessories Web Site

www.impact-test.com

User Guide

User Guide

User Guide

DO NOT ATTEMPT TO OPERATE THIS EQUIPMENT BEFORE READING AND COMPREHENDING THE MANUAL IN ALL ITS PARTS

USERS

- MACHINE MANUFACTURERS
- DRAUGHTSMEN
- OPERATORS
- MAINTENANCE WORKERS
- ANY OTHERS



WARNING

BEFORE USING THE ASPIRATOR SUPPLIED WITH THE SIEVE SHAKER, CHECK THAT THE PAPER AIR FILTER HAS BEEN TAKEN OUT FROM THE ASPIRATOR; IF IT IS STILL PRESENT, REMOVE IT.

BEFORE USING THE APPLIANCE, READ THE CHAPTER "SUGGESTIONS FOR THE ASPIRATOR USAGE" AND CONSULT THE OPERATING INSTRUCTIONS OF THE ASPIRATOR.

CONTENTS

Chapter 1	GENERAL INFORMATION
1.01	Warnings
1.02	WARNING and DANGER indications - SIGNS
1.03	Aim of the Instructions Manual
1.04	Modifications and enclosures of the Instructions Manual
1.05	N/A
1.06	EC Stamp
1.07	Usage
1.08	Operators
1.09	Storage
1.10	Transportation and Movement
1.11	Packaging removal
Chapter 2	TECHNICAL CHARACTERISTICS
2.01	General machine description
2.02	Dimension and weight
2.03	Electrical supply
2.04	Noise
Chapter 3	GENERAL SAFETY NORMS
3.01	General norms
3.02	Machine Safety devices and protection
3.03	Dangerous parts and residual risks
Chapter 4	INSTALLATION INSTRUCTIONS
4.01	Location
4.02	Electrical connection
Chapter 5	IN FUNCTION - USE
5.01	Set up
5.02	How to connect the aspirator to the sieve- shaker
5.03	Machine calibration – Meters - Indicators
5.04	How to place the material to be sieved
5.05	Switching the equipment on
5.06	Control panel adjustment
5.07	Vacuum adjustment
5.08	Emergency stop
5.09	Start up after emergency
5.10	Switching off
5.11	Test Start Up
5.12	Suggestion for the aspirator usage
5.13	Practical example of usage
Chapter 6	MAINTENANCE
6.01	Ordinary Maintenance
6.02	Extraordinary Maintenance
6.03	Authorised maintenance centres
Chapter 7	GUIDE TO RECOGNISING DAMAGE AND ANOMALIES
Chapter 8	SPARE PARTS
Chapter 9	INACTIVITY
Chapter 10	DECOMMISSIONING THE MACHINE

Chapter 1 GENERAL INFORMATION

1.01 WARNINGS

The manufacturer does not accept any responsibility for direct or indirect damage to people, things or animals and use of the appliance in different conditions from those foreseen.

The manufacturer reserves the right to make changes to the documentary information or to the appliance without advance notice.

Check the machine responds to the standards in force in the state in which it has been installed.

All operations necessary for maintaining machine efficiency before and throughout use are the operator's responsibility.

Carefully read the entire manual before operating the machine.

It is vital to know the information and limitations contained in this manual for correct machine use by the operator.

Interventions are only permitted if the operator is accordingly competent and trained.

The operator must be knowledgeable about machine operations and mechanisms.

The purchaser must ensure that operators are trained and aware of all the information and clarifications in the supplied documentation.

Even with such certainty the operator or user must be informed and therefore aware of potential risks when operating the machine.

Safety, reliability and optimum performance is guaranteed when using original parts.

Any tampering or modifying of the appliance (electrical, mechanical or other) which has not been previously authorised in writing by the manufacturer is considered abusive and disclaims the constructor from any responsibility for any resulting damage.

All necessary operations to maintain the efficiency of the machine before and throughout use are the responsibility of the user.

1.02 WARNING AND DANGER INDICATIONS - SIGNS

The machine has been designed and constructed according to the current norms and consequently with mechanical and electrical safety devices designed to protect the operator or user from possible physical damage. Residual risks during use or in some intervention procedures on the device are however present. Such risks can be reduced by carefully following manual procedures, using the suggested individual protection devices and respecting the legal and safety norms in force.

This manual includes "Warning" and "Danger" indications in relevant chapters. These indications are shown with the words "Danger" or "Warning" in bold font and uppercase to make them highly visible.

"**WARNING**" indicates that machine damage could be caused should indications be ignored.

"**DANGER**" indicates that machine damage and/or injury to the worker could be caused should indications be ignored.

"**DANGEROUS ZONE**" indicates any zone inside or in the proximity of the appliance in which a person is exposed to the risk of injury or damage to health.

1.03 AIM OF THE INSTRUCTIONS MANUAL

This manual has been edited with the aim of providing all machine operators with all the necessary information on installation, use and maintenance from production to scrapping in as comprehensive and clear manner as possible.

All the procedures useful for any foreseeable emergency situations have been listed by the manufacturer and can be verified during use.

Operators, for whom this manual has been written, due to their competence must give instructions or operate the machine themselves.

The instructions manual must be carefully consulted by laboratory or site safety managers, equipment operators and any internal and external maintenance workers.

The manual is integral to the product and refers to this appliance only.

The manual must be safeguarded and always kept near the equipment so that it can be easily consulted whenever necessary.

IMPORTANT: The manual does not substitute the experience and technical training of the worker but must be considered a guide for carrying out its functions.

Furthermore all the norms and rules the operator should be aware of or consult for correct use of the machine and/or test performance can be found in the manual.

This responsibility is entrusted to the installer and Laboratory or Site Manager where the machine is installed.

The Constructor is available to provide further information.

1.04 MODIFICATIONS AND ENCLOSURES OF THE INSTRUCTIONS MANUAL


This manual reflects the state at the time of the launch of the machine on its market. If any modifications, improvements or adjustments have been made since machine supply the Manufacturer does not have to intervene on the marketed machine and will not consider the machine or the manual deficient or inadequate.

1.06 EC STAMP

SEE EC DECLARATION

1.07 USAGE

The air jet machine is used for sieving powder and dry grain products by obtaining sieving results between 5 to 4000 microns. Its working foundation is based on the use of air that tug thin particles to make them go through the sieve. This effect is made through a vacuum machine that provokes a controlled decrease of pressure.

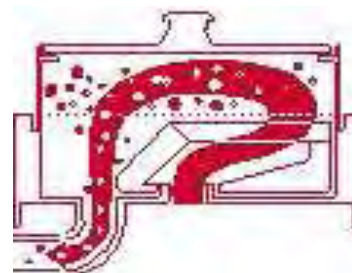

STANDARD EQUIPMENT

- Sieve shaker
- Supplying cable
- Plexiglass cover
- Little plastic hammer
- Instructions manual
- Vacuum cleaner and connection pipe

This appliance is for the exclusive use which it has been conceived for. Any other use is considered improper and therefore negligent. Machine use is allowed only in places free from danger of explosion or fire. During operation check for conditions of danger. immediately stop the machine should it be working irregularly, and consult the authorised dealer's Sales Service department. It is the Client's responsibility to verify at the time of installation and use that no conditions of use arise which are different to those indicated. Refer to the Constructor when in doubt.

1.08 OPERATORS

	DANGER WARNING	The use, transportation, installation, maintenance, demolition and disposal of the appliance are only permitted to "QUALIFIED PERSONNEL". This manual is exclusively aimed at "QUALIFIED PERSONNEL" and contains the necessary information for machine use.
--	-----------------------	---

"QUALIFIED PERSONNEL" means people who, due to their training, experience and education, as well as knowledge of the relevant standards, limitations and measures, have been authorised by the "PLANT SAFETY MANAGER" to carry out any necessary activity and are able to recognise and avoid any possible danger.

The manufacturer recommends that the instructions, procedures and recommendations in this manual and the work safety legislation in force be scrupulously adhered to, even with the use of appropriate protection devices (whether individual or part of the machine).

Knowledge and respect of the instructions, safety warnings and danger in this manual are all necessary for installation, operation, management and machine maintenance with a minimal risk.

The "PLANT SAFETY MANAGER" has the following responsibilities and duties:

- To know the machine functions, its commands, safety and protection devices, possible dangers of use and all the information in this manual in detail. This knowledge can only be gleaned from detailed reading of this manual.
- To know the safety legislation in force in detail in order to operate the machine

- To recognise the “QUALIFIED PERSONNEL” for transportation, handling, installation, use, maintenance, disposal, etc.
- Correctly train and educate the “QUALIFIED PERSONNEL” before allowing them access to the machine. The personnel must also be exhaustively trained with regards to the machine’s protection devices.
- Ensure the machine’s safety devices are not tampered with or removed and are checked on a daily basis. Provide the operator appropriate individual protection devices according to the laws in force.
- The constructor is available for clarification, assistance and training and declines all responsibility for damage to things or people resulting from improper, incorrect and negligent use by untrained personnel.

1.09 STORAGE

	WARNING	The appliance must be stored and conserved in the original packaging and in a closed environment, protected from atmospheric agents with a minimum temperature of -15C°, and a maximum of +60C° and a maximum humidity of 70 %.
--	----------------	---

1.10 TRANSPORTATION AND MOVEMENT

	WARNING	In order to avoid irreparable machine damage, move with care, do not overturn, protect from rain, do not stack, protect the packaging and its contents from bumps and sources of heat.
--	----------------	--

During transportation and movement it is important to avoid bumps, overloading with other packages, exposure to freezing or heating atmospheric agents, or any other potentially harmful condition to the device, things or people. Machine transportation and movement must be entrusted to Qualified Personnel who can ensure correct movement.

		DANGER WARNING	Do not transport or move the product should it be impossible to respect the conditions on the packaging or there be any doubts. Request information from the constructor.
--	--	-----------------------	---

1.11 PACKAGING REMOVAL

After removing the packaging check the machine is complete and that there are no visibly damaged parts. **DO NOT USE THE MACHINE** and refer to the constructor when in doubt.

	DANGER	The components used for packaging (plastic bags, polystyrene, nails, screws, wood, etc) must be kept out of reach of children, as they are sources of danger. These components should be placed in the appropriate containers.
--	---------------	--

	WARNING	In order to avoid bumps and overturn adopt the normal and logical precautions. Before disposing of the packaging check all machine components such as accessories, utensils, instructions, documents etc have been removed.
--	----------------	---

Chapter 2 TECHNICAL CHARACTERISTICS

2.01 GENERAL MACHINE DESCRIPTION

LID A1

It is used to seal the sieve **A2** and the material that have to be sieved during the test in order to avoid material's outflow.

SIEVE A2 (it's an accessory not supplied with the instrument).

GENERAL SWITCH A3

The switch enables or disables all the sieving-shaker functions, including the aspirator functioning. Next to the general switch there is the supplying voltage tap with the appropriate connection cable supplied with the equipment.

ELECTRIC TAP FOR ASPIRATOR A4


Tap to which is connected the electric supplying of the aspirator.

MOUTH PIECE OF THE ASPIRATOR PIPE A5

The aspirator pipe is connected to this mouth piece; the aspirator is used to create the vacuum

REGULATION COUPLING OF THE VACUUM A6

After starting the aspirator, it is possible to increase or decrease the generated vacuum degree by screwing and unscrewing this regulator device.

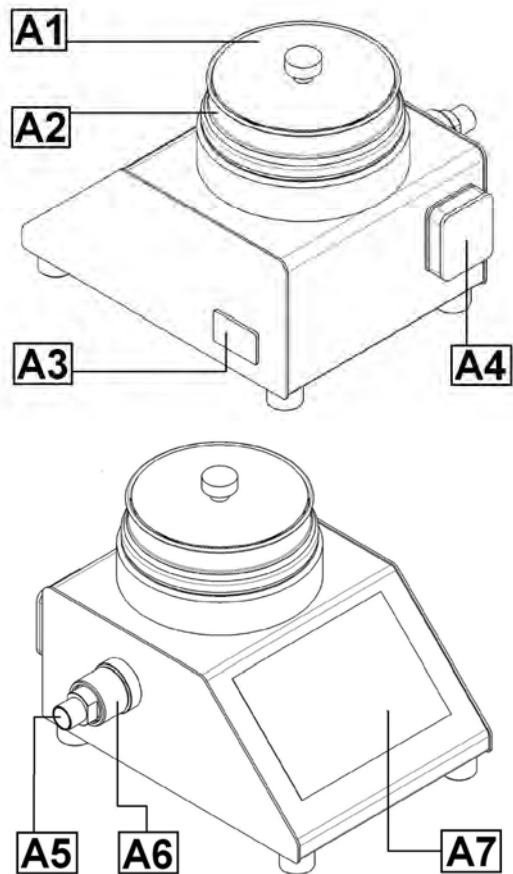

2.02 DIMENSION AND WEIGHT

LENGTH	430 mm
WIDTH	407 mm
HEIGHT	340 mm
WEIGHT	20kg circa

2.03 ELECTRICAL SUPPLY

VOLTAGE	230 V
PHASE NUMBERS	1
FREQUENCY	50Hz
SIEVING MACHINE POWER	100 WATT
ASPIRATOR POWER	1300W

2.04 NOISE

The air noise emission levels shown do not necessarily imply the levels of exposure to the worker.

The levels of exposure to the operator are obviously linked to the emission levels of the appliance; however other factors influence the levels of exposure to the operator: length of exposure, environmental characteristics, the presence of other machines etc.

The appliance emission levels allow anyway an estimate to be carried out on the dangers due to noise.

**DANGER**

Continuous use of the appliance and machines predictably present in the installation environment could cause a heightened daily personal exposure to noise.

If the daily personal exposure is equal to or more than 85 dB (A) it is advisable to use the Individual Protection Devices (protective headphones, plugs, etc.). If the daily personal exposure is equal to or more than 90 dB (A) it is compulsory to use Individual Protection Devices (protective headphones, plugs, etc.). For further information consult the standards in force in the country of installation.

Average continuous equivalent acoustic pressure level L_{Aeq} to the distance 1.5 m 61 (A):
Level of acoustic power emitted by the LWA device xx dB (A):
Standard EN ISO 3746

Chapter 3 GENERAL SAFETY STANDARDS**3.01 GENERAL STANDARDS**

To ensure the safety of machine operators:

- Any tampering with the appliance not pre-emptively authorised by the manufacturer exempts the manufacturer from any responsibility for damage caused by or to it.
- The removal or tampering with safety devices entails a violation of the safety standards.
- Machine use is only allowed in areas where there is no risk of explosions or fires.
- Only the original fittings can be used. The use of unoriginal fittings exonerates the manufacturer from all responsibility.
- Check the appliance is in ideal working conditions and that its parts are not worn or faulty before Carry out all necessary maintenance
- Do not wear loose clothing, ties, chains or anything else which could become caught in the frame or other moving parts of the appliance.
- Be aware of the danger of electrical shocks from direct or indirect contact due to unforeseen electrical faults.
- Do not subject the appliance to violent impact.
- Do not expose the appliance to fire, welding sparks or extreme temperatures.
- Do not bring the appliance into contact with corrosive substances.
- Do not wash the appliance with jets of water.
- Check the workspace around the machine is clear from potentially dangerous objects.
- The machine operator must wear appropriate work clothing such as protective glasses, gloves and mask in order to avoid damage from, for example, harmful dust projection. Wear a lower back support when lifting heavy parts. There should be no hanging objects such as bracelets or otherwise, long hair should be protected with relevant precautions, shoes must be appropriate for the type of operation to be carried out.

DURING USE

When operating check there are no conditions of danger. Immediately stop the machine when it is functioning irregularly. Contact the authorised Sales Service department.

- For the operator's safety do not touch any part of the appliance when testing and use the appropriate individual protection devices in order to keep the operator safe.

3.02 MACHINE SAFETY DEVICES AND PROTECTION

DEFINITION: Protections are all the safety measures that consist of the use of specific technical means (repairs, safety devices) to protect people from dangers which cannot be limited reasonably in design.

**DANGER**

Tampering with the protections or any appliance modification could cause risks to users or other exposed people.
The manufacturer does not assume any responsibility for direct or in direct damage to people, things or animals following tampering with the protections.

The sieving machine is provided with a general switch **A3**, that works also as emergency switch.

3.03 DANGEROUS PARTS AND RESIDUAL RISK

During some procedures the operator could face some risks of danger.
The risks can be eliminated following carefully the procedures written in this manual and using suitable safety devices

RISK OR DANGER	PROTECTION DEVICE
FINGER SQUEEZE	REINFORCED GLOVES
LOSS OF GRANULATED MATERIAL	GLASSES

DURING THE USE

In order to assure the operator safety, it is recommended to:

1. Do not remove the lid and the sieve during the functioning of the sieving machine
2. Do not ever switch the sieving machine on if the lid and the sieve have not been positioned
3. Carry out the equipping and maintenance operations only when the feeding cable is disconnected from its own supplying tap.

Chapter 4 INSTALLATION INSTRUCTIONS

4.01 LOCATION


The equipment must be placed in an ideal position and environment for the use it has been conceived for (laboratory use and protected from atmospheric agents) and that the machine is placed by a qualified operator.

ALLOWED TEMPERATURE:	from +5°C to +40°C
ALLOWED RELATIVE HUMIDITY:	from 30% to 70%
MAXIMUM HEIGHT OVER SEA LEVEL:	1000 m


GENERAL ADVICE

- The machine must be installed in an area which allows ease of access to all parts so that maintenance may be carried out.
- Unauthorised people and objects which could be potential sources of danger must not be permitted in the area surrounding the machine.
- Do not position the equipment near instruments or appliances which could produce vibrations.


The appliance rests on rubber feet and had to be placed on a plane surface for a better stability, otherwise it may capsize and cause injuries.

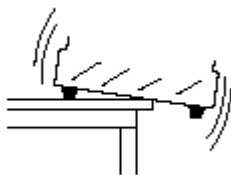

For an easy use it is advisable to place the appliance at about 700 mm height from the ground surface.
The required minimum space for the installation is 430x407x340. The control panel preferably has to be placed and fixed at such a distance that doesn't interfere with the equipping and sample loading operations.

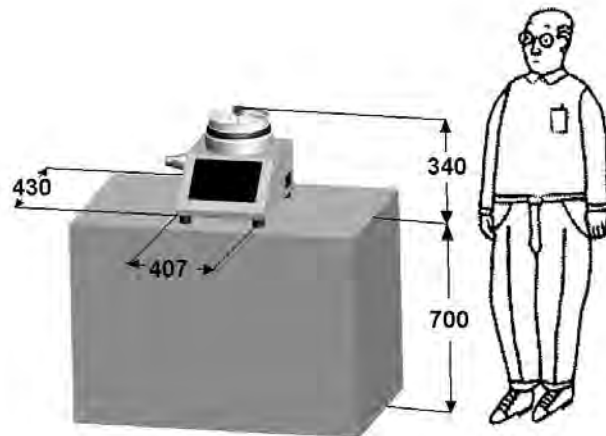

4.02 ELECTRICAL CONNECTION


	<p>DANGER</p>	<p>Wiring of the electrical system must be carried out by qualified personnel. Before wiring consult the electric plan linked to the instructions manual and the registration plate on the machine for information regarding supply, frequency and nominal current.</p> <p>Connect the earthing system via the PE terminal (yellow-green) before any other connection.</p> <p>Apply a knife switch at the top of the connecting cable of the machine to the power system.</p> <p>The knife switch must be combined with a safety device against the overload with a differential switch (safety switch).</p> <p>The technical features of the safety device must be in accordance with the standards in force in the country where the machine has been installed.</p>
---	----------------------	--

ELECTRIC TOLERANCES:

- Real voltage $\pm 10\%$ of the nominal one
- Frequency: $\pm 1\%$ of the nominal one in a continuous way
 $\pm 2\%$ of the nominal one for a short period
- The harmonic distortion of the sum from the second to the fifth harmonics not more than 10% of the total voltage as a real value between the conductors. A further distortion of 2% is accepted for the sum from the sixth to the thirtieth harmonics of the real total value between the conductors.
- With reference to the voltage imbalance of the three-phase voltage, the inverted sequence component and the zero sequence component must not be more than 2% of the direct sequence component of the voltage.
- The voltage pulses must not last more than 1,5 ms with an up/down time between 500 ms and 500 ms and a peak value not higher than 200% of the real value of the nominal tension.
- The electric supply must not be interrupted or zeroed for more than 3 ms at any time. Between two interruptions it must not take more than 1 s.
- The interruptions must not overcome 20% of the tension peak for more than one cycle. Between two interruptions it must not take more than 1 s.



The manufacturer assumes no liability for any damages to people, things and animals caused by the non-compliance of the above instructions

Chapter 5 IN FUNCTION - USE

	<p>DANGER WARNING</p>	<p>Before setting the machine in motion it is essential that the Operator and Safety Manager have read the Instructions Manual and understood all parts of the machine and activities linked to it (Risks, Dangers, Functionality, Operation, Protections, Commands, etc.)</p>
---	------------------------------	--


5.01 SET UP

Before using the appliance, check that the distance between the sieve bottom and the aspirator mouth is between 2 and 4 mm; this distance has been already checked in the factory during the production phase.

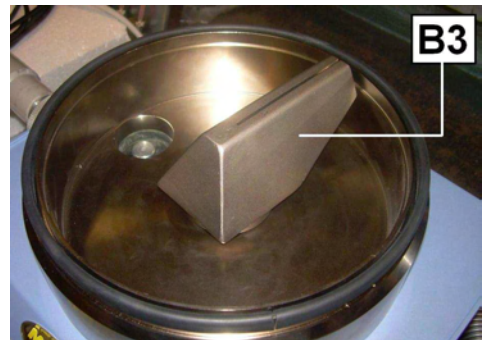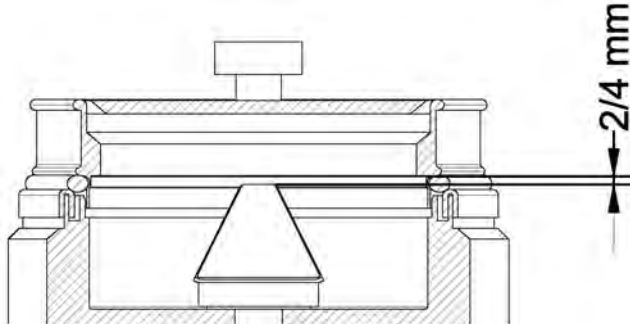

It is suggested to make a periodical inspection, especially if several types of sieves are used, by following these instructions:


1. Unscrew the aspirator mouth **B3**
 2. The rod **B2** is supplied with a regulation screw **B1** on its end **that determines the aspirator mouth's height compared to the sieve.** By tightening the screw the distance between the sieve and the aspirator's mouth increases, while by unscrewing it the distance decreases.
- ATTENTION: the opening of the aspirator's mouth have to be free of deposits.

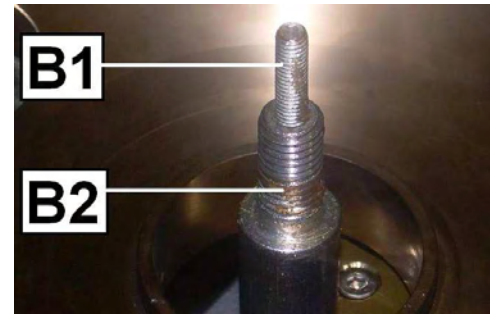

5.02 HOW TO CONNECT THE ASPIRATOR TO THE SIEVE- SHAKER

In order to start the sieving it is necessary to connect the aspirator to the sieve shaker as explained in the following instructions:

- connect the aspirator's pipe to the mouth piece **A5** of the sieve shaker (pic.3).
- connect the aspirator's feeding cable to the input **A4** that is placed on the back side of the sieve shaker.

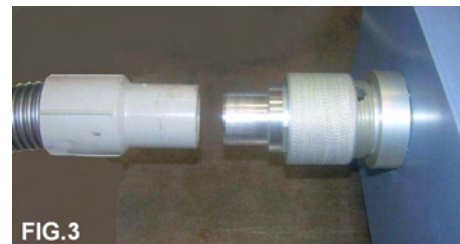

5.03 MACHINE CALIBRATION – METERS - INDICATORS

The machine is checked in the factory, using sample equipment periodically checked by officially recognised institutes. These checks cannot guarantee that the machine, meters and indicators will provide accurate values and results conforming to the standards in force in the countries the machine has been installed and used in. Normally such norms envisage calibration check after every movement. In order to obtain correct values and results it is therefore VITAL that the operator, once the machine has been installed and set up and before official tests, has an officially recognised body check the machine characteristics, its calibration and results/values reliability. The manufacturer is exempt from all responsibility in the case of direct and indirect damage from use of the machine without official approval by the relevant bodies.

As explained before, the sieve shaker is supplied already calibrated; in case it was necessary or it was required to check that the visualized vacuum value is correct, you have to contact the technical assistance service.


5.04 HOW TO PLACE THE MATERIAL TO BE SIEVED

The material which have to be sieved must be put in the upper sieve; after that, cover the sieve with the lid **A1**. If there is an high quantity of material, the sieving will take longer and the test will be more accurate.

The maximum material size must be checked. The material (that must be granulated) have to move on the sieve when the aspirator mouth is working.

The weight have to be in compliance with the kind of test to be executed. Our general recommendations for Ø 200 mm are:

- Normal condition. Sieves with a 0,040 mm opening: 20 gr.
- Heavy materials with good smoothness: up to 50 gr.
- Light materials hard to be sieved and sieves with opening less than 0,040 mm: 10gr.

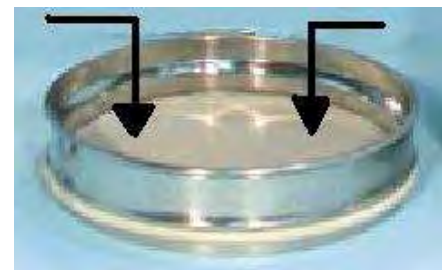

5.05 SWITCHING THE EQUIPMENT ON

To switch the sieve shaker on press the general switch **A3** and place it on the "I" position.

5.06 CONTROL PANEL ADJUSTMENT


The control panel is composed by 4 keys and one luminous display; it allows to select and visualize the main functioning parameters of the sieve shaker.

CONTROL PANEL OPERATING FEATURES

The control panel is composed by:

- **A display C1** composed by 4 figures, divided in 2 groups of 2 figures each; the 2 figures on the left shows the reached vacuum value (expressed in millibar), the 2 figures on the right shows the selected total time of the sieve-shaking.
- **A starting key C2**: by pressing it the sieving starts working.
- **A stop key C3**: it is used to stop the timer and sieving phase.
- **2 keys C4**: used to increase and decrease the sieving time

Every time that a key is pressed a bip will be given out.

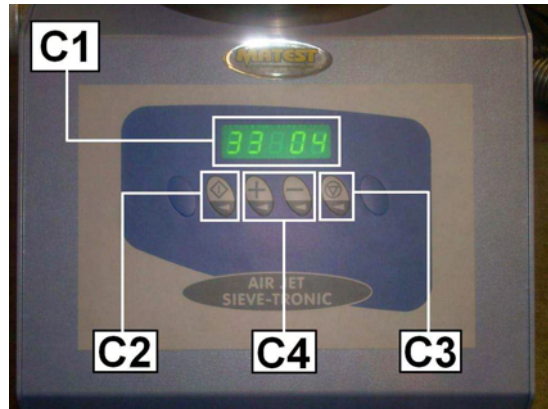

TIMER'S ADJUSTEMENT

After switching the sieve-shaker on and before beginning the sieving by pressing the 2 keys **C4**, it is possible to increase and decrease the sieving time. After having selected the desired time, press **C2 once** to save the selected value.

ATTENTION: it is necessary to press C2 once to save the setting every time the sieving time has been modified.

START-UP

Press **C2** to start-up the sieve-shaking; the timer will automatically begin the count-down of the selected time and will stop at the end; when the sieve-shaking ends the previously selected time will automatically appear on the screen.

NORMAL STOP

It is possible to stop the sieve-shaking before the end of the selected time, by pressing **C3**; **when the sieve-shaking is restarted, the previously selected time will automatically appear on the screen**

HOW TO STOP MOMENTARILY THE SIEVE SHAKING (PAUSE FUNCTION)

During the sieve-shaking it is possible to suspend it momentarily by pressing **C2** once: that enables the operator to run settlement operations of the material as well as equipping operations; the pause is pointed out by the flash of the figures on the display, showing the missing time.

To re-start the sieving press **C2** again; **in this case the count-down will start again from the point in which it had been previously stopped.**

HOW TO RE-SET UP THE TIMER DURING THE SIEVE-SHAKING

Press **C2** During the test to modify the remaining sieving time; the sieving will be immediately stopped; the suspension is pointed out by the flashing of the figures indicating the remaining time; at this point it is possible to adjust the timer using the 2 keys **C4**.

After having reselected the sieving time, press 2 times **C2** to restart the sieve shaking (**once to save the selected time and once to re-start the test**).

5.07 VACUUM ADJUSTMENT

To adjust the vacuum produced by the sieve-shaker, it is necessary to roll **A6**; unscrew the coupling to reduce the depression and screw it to increase the depression.

The increase and the decrease of the vacuum produced by the aspirator will be displayed on the screen.

In accordance to UNE-EN 933-10 Standard "Test for the determination of arid geometric properties" the recommended pressure is $3,0 \pm 0,5$ kPa (-30 millibar). Anyway we suggest to run more tests in order to be able to determine the correct pressure for different kinds of materials. The too soft materials (easy to be pulverized) must be sieved with the possible lowest pressure.

5.08 EMERGENCY STOP

In case of emergency it is possible to stop the sieve-shaker by pushing the general switch **A3** and by placing it on 0 position.

5.09 START UP AFTER EMERGENCY



**WARNING
DANGER**

Before re-starting the appliance determine and eliminate the causes of the emergency stop

To re-start the sieve-shaker push the general switch **A3** and place it on "I".

5.10 SWITCHING OFF

Press **C3** to stop the sieving machine

5.11 TEST START UP

Before using the appliance regularly check it is working correctly by carrying out at least one complete empty cycle according to the previous instructions.

Should there be any problem consult the chapter "DIAGNOSTICS".

If the instructions in this manual do not provide the solution to the problem, contact Sales Assistance.

5.12 SUGGESTIONS FOR THE ASPIRATOR USAGE

Some suggestions are listed as follows; they allow the operator to use the aspirator correctly:

- Remove or do not assemble the paper filter on the aspirator in order to use it correctly, but leave the plastic bag.
- During the test, every time that the vacuum created by the aspirator is not enough to satisfy the requirements declared by the Standard, activate the control of the filter cartridge automatic cleaning (the aspirator's filter cartridge could be dirty); in this way the aspirator will work at its maximum potentialities.
- If the vacuum created by the aspirator will not be enough even after the filter cleaning operation, it is necessary to replace the filter cartridge placed in the sieve shaker. Consult the instructions enclosed to the aspirator in order to replace the cartridge filter of the sieve shaker.



WARNING

NEVER SUCK IN WITHOUT THE CARTRIDGE FILTER

5.13 PRACTICAL EXAMPLE OF USAGE

Hereunder we describe a "standard-procedure" which allows even to an operator without a wide experience to carry out a test.

Naturally the operator experience will allow to optimise the machine performances on the basis of the customer needs

In order to have several points in the grinding material curve, start the test with the sieve with smaller openings and sieve the material that has not been passed through the sieve with larger openings. Follow the same procedure using other openings and repeat it depending on how many divisions the operator would like to obtain.

The hygroscopic products absorb the humidity of the air during the sieving procedure. This could be the reason of some measuring mistakes.

To avoid this kind of mistakes it is suggested to dry the material before and after every test.



DANGER

Do not attempt to use this machine without reading and fully understanding this manual in all parts.

**WARNING**

Do not assemble the paper filter but only the plastic bag for a correct use of the aspirator.

1. Check the distance between the sieve and the aspirator mouth as described in chapter “**SET UP**”
2. Position the material to be sieved as described in chapter “**HOW TO PLACE THE MATERIAL TO BE SIEVED**”
3. Connect the aspirator to the sieve shaker as explained in chapter “**HOW TO CONNECT THE ASPIRATOR TO THE SIEVE- SHAKER**”
4. Switch the sieve-shaker on by pressing the general switch **A3**.
5. Adjust the timer as described in chapter “**TIMER’S ADJUSTMENT**”. In accordance to the UNE-EN 933-10 Standard, 2001 edition, the sieve shaking time can be considered concluded when the weight does not change more than 0,1% in one minute.
6. Start the instrument by pressing **C2**
7. During the sieve-shaking process it is advisable to hit the cover with the plastic hammer to remove the dust of the material. It is possible to stop the machine under vacuum and the air sieve-shaker to remove with a brush the residual materials that may have produced.
8. Wait or the end of the sieve-shaking time selected on the timer, when the time is expired the sieving will stop (to stop the test before the selected time press **C3**).
9. When the sieving ends, it is necessary to weigh the material that is still on the sieve. To avoid loss of material, brush the inner surface of the cover and brush accurately the openings when the pulverized granules are put on another sieve.

Chapter 6 MAINTENANCE**6.01 ORDINARY MAINTENANCE****WARNING
DANGER**

Do not perform maintenance – interventions on the machine which have not been quoted and described in this instructions manual without first contacting the manufacturer.
Periodically clean all machine parts and oil the unpainted parts in order to preserve the machine and its efficiency.
Avoid the use of solvents which damage paint and parts in synthetic material.

How to clean the sieve-shaker:

- To remove the aspirator’s mouth, pull it upwards.
- Connect the sieve-shaker and clean the aspirator’s mouth with a brush and the sieve- shaker’s body and its inner part with a vacuum cleaner.

ATTENTION: do not turn the aspirator’s mouth **B3** with hands

How to clean the aspirator:

Consult the chapter “**SUGGESTIONS FOR THE ASPIRATOR USAGE**” for the aspirator cleaning operations.

How to clean the sieves:

Clean carefully the lower part of the sieve with a soft brush. Do not push and do not hit the sieve’s mesh.
The most efficient way to clean the sieves is an ultrasonic cleaner. An ultrasonic cleaner is the fittest for this use. To increase the effectiveness of the cleaner, use a water detergent.

6.02 EXTRAORDINARY MAINTENANCE

For extraordinary maintenance operations refer directly to the Manufacturer.

6.03 AUTHORISED MAINTENANCE CENTRES



For information on the nearest authorised help centre it is essential to contact the manufacturer.

Chapter 7	GUIDE TO RECOGNISING DAMAGE AND ANOMALIES
------------------	--

This chapter presents and discusses all the simple problems which could occur during machine use. The appropriately qualified, professional personnel must carry out all the maintenance procedures, check and control, as well as all the repair operations on parts of the machine or the electrical system. Contact Technical Sales Assistance for any other problem not listed on the previous table or should the malfunctioning persist after the intervention of the operator in accordance with the previously mentioned courses of action.

PROBLEM	POSSIBLE CAUSE	REMEDY
When pressing A3 the control panel does not switch on	Fuses are broken	Replace them
	Lack of supply	Check the connection to the power grid
When pressing A3 the control panel switches on but the aspirator's mouth does not turn.	Damage in the electric system	Check the electric system
	Ratiomotor malfunctioning	Ratiomotor replacement
It is not possible to reach the vacuum value as required by the Standard	Damage of the vacuum sensor	Replace the sensor
	Blow-by from the suction equipment	Find the 3 filaments and remove them
	The aspirator does not work properly	Clean the filters; look at the "SUGGESTIONS FOR THE ASPIRATOR USAGE" chapter.

Chapter 8	SPARE PARTS
------------------	--------------------

 	WARNING DANGER	<p>Only original spare parts can be used. Use of unoriginal spare parts exempts the manufacturer from all responsibility. Procedures for substitution of spare parts will be provided by the manufacturer along with the part. For spare parts contact the manufacturer's Sales Service department.</p>
---	---------------------------	---

SIEVE A058-12:

Sieve Ø200mm, stainless steel frame and nylon cloth , openings from 5 to 71 microns (when ordering please specify cloth opening).

SIEVE A058-13:

Sieve Ø200mm, stainless steel frame and cloth, openings from 75 to 4000 microns (when ordering please specify cloth opening).

FILTER CARTRIDGE A058-14:

Spare part for the aspirator filter cartridge replacement.

PLASTIC BAGS A058-15:

Spare part : pack of 5 pieces.

Chapter 9	INACTIVITY
------------------	-------------------

Ensure all machine parts are in safe working order before operating it again should the machine be inactive for a long period of time. When in doubt contact the Manufacturer

Chapter 10 DECOMMISSIONING THE MACHINE

Should it be decided that the machine is to be no longer used, proceed as follows:

- Disconnect the electrical supply network by removing the connecting cable therefore making it unusable.
- Make the potential sources of danger harmless, such as sharp or protruding parts.
- Dismantle the machine; divide it into similar parts and dispose of according to the standards in force.

Recycling notice for the disposal of electrical and electronic devices



This symbol, shown on the device or on the package and/or the documentation, suggests that the device should not be disposed together with other home garbage at the end of its life cycle.

To avoid further environment, or health-care damage, caused by the unsuitable disposal of garbage, the user should separate this device from other different types of garbage and recycle it in responsibly to avoid the reuse of material resources. Users must take care at the disposal of the equipment by taking it to the nearest recycling site for appropriate recycling treatment for electrical and electronic devices. Gathering and Recycling deplete devices allow the preservation of natural resources and grant them the adequate treatment by respecting health and environment.

For further information on your local recycling site please contact your local council or city waste treatment department. The developer, as producer of electrical and electronic devices, will provide to finance the recycling and treatment services for deplete devices that will come back through these recycling sites, according to the local statement.