

Tabletop Aluminum Cap Sealer - Manual

BAL-5020 BAL-5015 BAL-5013

Dimension

Height : 35cm

Weight : 5 kg


Size of clip : BAL-5020 20mm
BAL-5015 15mm
BAL-5013 13mm

Introduction:

Portable and convenient laptop sealing machine specialized in aluminum lips. Competitive price, easy operating, can achieve safety speed and time saving requirements.

Application:

Injection pharmacy, medicine, cosmetic, contact lenses...etc.

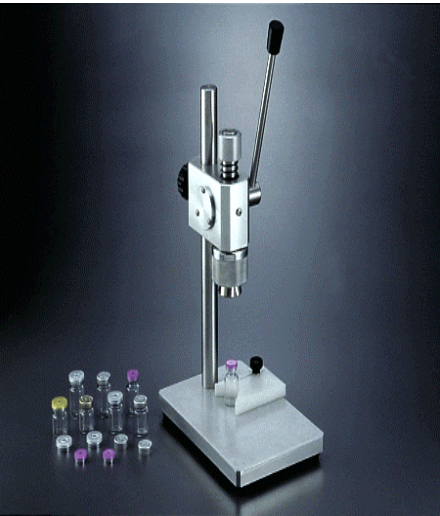

Accessories :

item	Name of article	No.	Quantity	Description
1	Manual		1 copy	
2	Special screw (round head)		3 PCS	backup
3	Special screw (flat head)		3 PCS	backup
4	Cross-screwdriver		1 PC	
5	Nut fixing tool		1 PC	
6	Accessory		1 tin	

Filler Co., Ltd.


Tel : (02)2593-1616 / 2591-6873

Fax : (02)2586-4587

www.filler.com.tw

E-mail: filler@filler.com.tw


Structure :

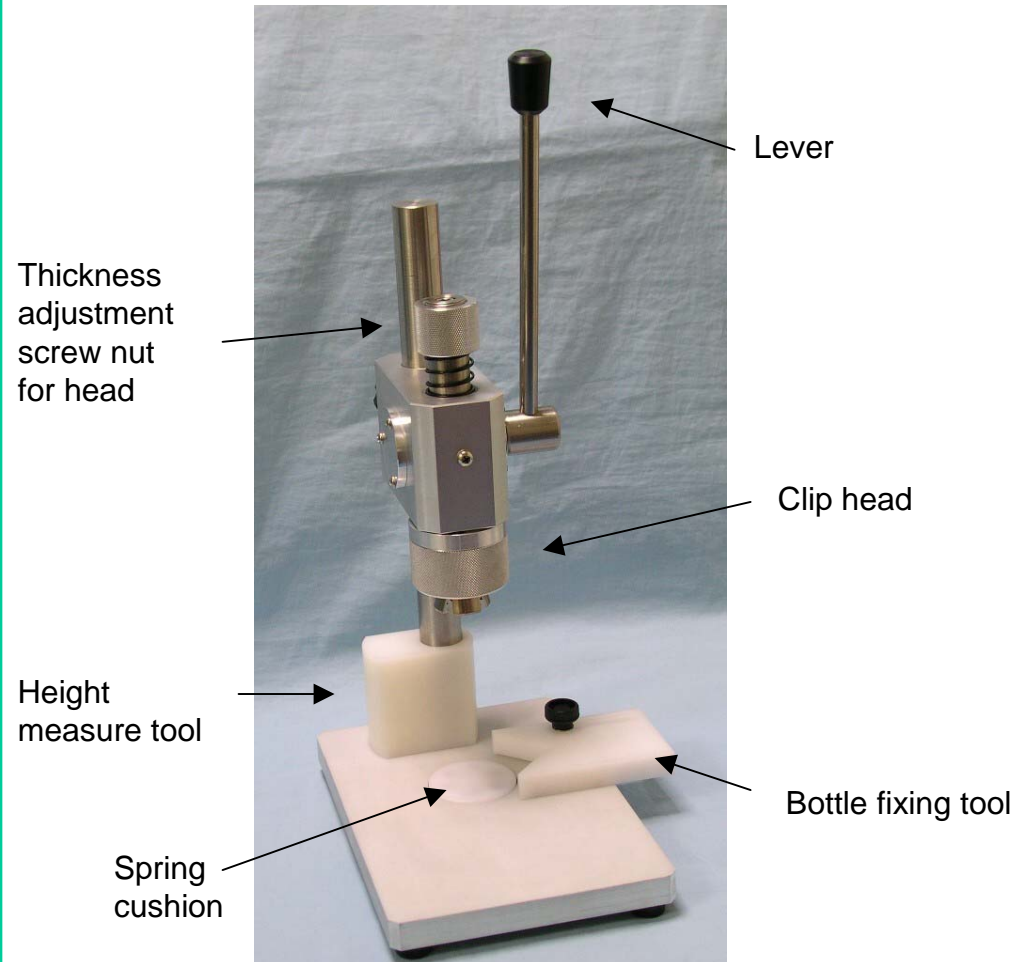


Note : Only 2 pieces of clip and above attach the tool and accessories.

Assembling Procedure :

1. First, screw the clip head(1) counterclockwise. Insert the nut fixing tool(2) for assistant.

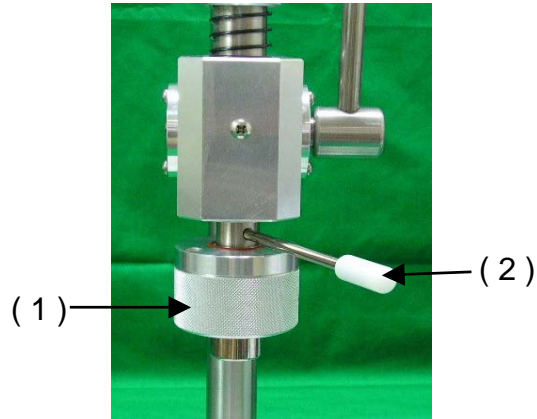

2. Press the lever(1) then screw (2) under the clip head by cross-screwdriver(3), assembling procedure completed.

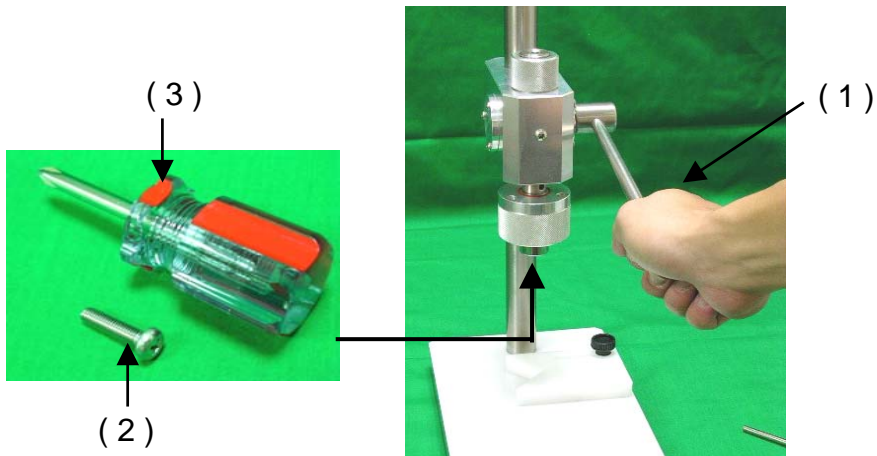

When the clip head can't suit the size of aluminum lip in normal use, adjust the nut(1) to loosen or fasten.

Screw the nut clockwise--loosen




Screw the nut counterclockwise--fasten




Disassembling Procedure : reverse step according to the assembling procedure.

Height Adjustment:

1. Put the bottle without the plug and cap under the height measure tool (1).
Loosen the screw (2) to put down the main machine on the height tool (3) then screw it on.

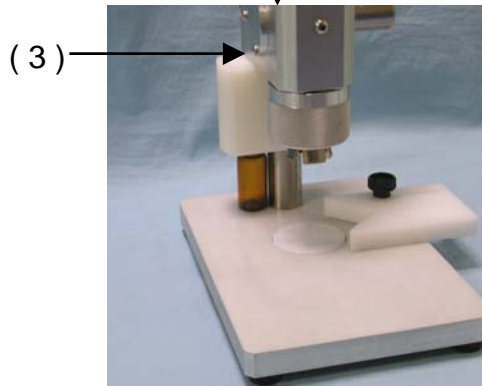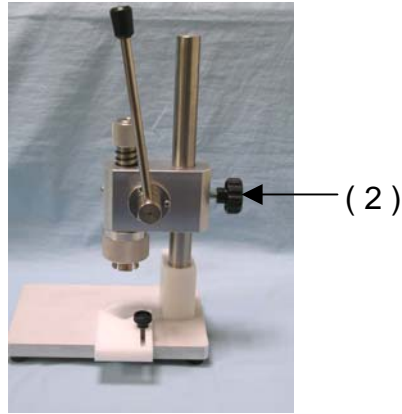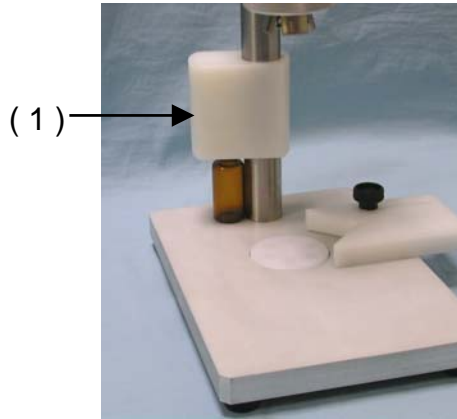

2. Put the bottle with the plug and cap under the clip head (1), please note the position should be in the center of cushion. Press the lever to seal the cap(2) then move the fixing bottle tool (3) to fix the bottle. Start the sealing process.

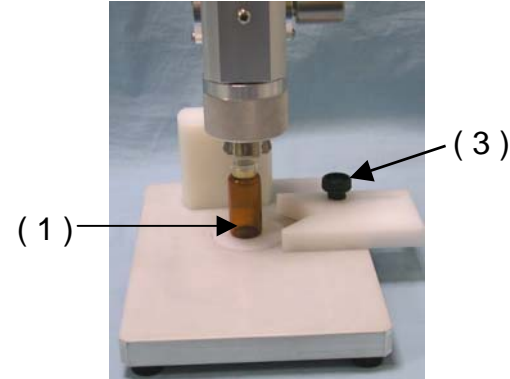

One (1) Year Limited Warranty - Worldwide

Filler Co. warrants their standard products against defects in their operation and non-consumable materials under normal use for a period of ONE (1) YEAR from the date of purchase by the original purchaser ("Warranty Period"). When a non-consumable hardware defect arises within the Warranty Period, a claim written in detail is to be received by Filler Co. within 30 days of the date the defect was discovered. At its option, Filler Co. will either (1) repair the hardware defect at no charge, using new or refurbished replacement parts, or (2) exchange the product with a product that is new or which has been manufactured from new or serviceable used parts and is at least functionally equivalent to the original product. This Limited Warranty does not apply to the damage caused by failure to either (1) follow instruction relating to the product's use, or (2) handle, fill, transmit, wash, clamp, and/or manufacture the materials that are feasible and agreed upon through product testing beforehand and under general industrial acceptance.

Except as provided in this Warranty and to the extent permitted by law, Filler Co. shall not be liable for any direct, special, incidental, or consequential damages of any kind, including, but not limited to any expenses for removal or reinstallation resulting from a defect; loss of use; loss of revenue; loss of actual or anticipated profits (including loss of profits on contracts); loss of the use of money; loss of anticipated savings; loss of business; loss of opportunity; loss of goodwill; loss of reputation; loss of, damage to or corruption of material; or any indirect or consequential loss or damage howsoever caused including the replacement of equipment and property, any costs of recovering, programming, or reproducing any material stored, filled, transmitted, handled, washed, clamped or used with Filler Co.'s products.

In practical terms:

- 1) We warrant this machine with quality of performing steady operation in normal use (based on existing industrial standard) according to its design, for one year counting from the date of purchase.*
 - 2) Parts that are considered consumables (such as Filling Pipe, Hand-held extension Pipe, Intake Pipe, etc.) are not covered under this warranty.*
 - 3) Parts & labor will be charged additional if the breakdown of machine is due to following:*
 - Unauthorized disassembling of the machine.*
 - Unauthorized operation based on instructions NOT included in User's Manual.*
 - Inappropriate power supply voltage applied.*
 - Damage from rats, insects, poor working condition, and natural disasters.*
 - Accident caused by failure in following instructions on assembling, disassembling, and cleaning procedures. (for example, sinking/soaking machine into water)*
-

* Proof of Purchase *

Owner's Name			
Address			
Article			
Purchase Date			
Dealer			Signature
Model			
TEL			
FAX			
Manufacturer	Filler Co.,LDT.		Signature
Model	BAL-5020		
TEL	886-2-2593-1616	886-2-2591-6873	
FAX	886-2-2586-4587		

1. We warrant the quality of this machine, in normal use, for one year from the date of purchase.
Consumables are not under warranty.
 2. Under warranty, partial material costs and service charges are necessary to repair the breakdown owing to :
 - a. Irregular use against the User's Manual
 - b. Unauthorized disassembling
 - c. Inappropriate power supply voltage applied
 - d. Private modification
 - e. Disasters, rats and insects
 - f. Personal factors (sinking / soaking into water) to cause severe damage
-