

Applied Biosystems S0LiD[™] 4 System

SOLiD[™] Bisulfite-Converted Fragment Library Preparation Protocol

> Bisulfite-Converted Library Preparation

Templated Bead Preparation Instrument Operation



For Research Use Only. Not intended for any animal or human therapeutic or diagnostic use.

This protocol is the proprietary material of Applied Biosystems, LLC or its affiliates and is protected by laws of copyright. The customer of the SOLiD™ System is hereby granted limited, non-exclusive rights to use this protocol solely for the purpose of operating the SOLiD™ System. Unauthorized copying, renting, modifying, or creating derivatives of this protocol is prohibited.

Information in this document is subject to change without notice.

APPLIED BIOSYSTEMS DISCLAIMS ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT OF INTELLECTUAL PROPERTY TO THE FULLEST EXTENT ALLOWED BY LAW, IN NO EVENT SHALL APPLIED BIOSYSTEMS BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF, WHETHER OR NOT FORESEEABLE AND WHETHER OR NOT APPLIED BIOSYSTEMS IS ADVISED OF THE POSSIBILITY OF SUCH DAMAGES.

TRADEMARKS:

The trademarks mentioned herein are the property of Life Technologies Corporation or their respective owners.

NanoDrop is a trademark of NanoDrop Technologies, LLC.

Agentcourt and AMPure are trademarks of Beckman Coulter, Inc.

Eppendorf is a trademark of Eppendorf AG.

Covaris is a trademark of Covaris, Inc.

TaqMan is a trademark of Roche Molecular Systems, Inc.

© Copyright 2010, Life Technologies Corporation. All rights reserved.

Part no. MAN0003271

Rev. date 2 December 2010

Table of Contents

	Required Materials	4
	Description of the Protocol	7
Metho	ods	9
	Before Starting	9
	Step 1. Shear the DNA	10
	Optional: Spike-In Lambda DNA Digested with Alu I	11
	Step 2. Prepare the Special dNTP Mixes	12
	Step 3. Prepare the Methyl-P1 Adaptor	13
	Step 4. End-Repair the DNA	15
	Step 5. Purify the End-Repaired DNA	16
	Step 6. Ligate Methyl-P1 and P2 Adaptors to the DNA	17
	Step 7. Purify the Adaptor-Ligated DNA	18
	Step 8. Nick-Translate the DNA	19
	Step 9. Purify the Nick-Translated DNA	20
	Step 10. Bisulfite Convert the DNA	21
	Step 11. Desulfonate and Purify the Converted DNA	23
	Step 12. Amplify the DNA	25
	Step 13. Purify the Amplified DNA	27
	Step 14. Size-Select the DNA	28
	Step 15. Quantitate the Library	30
	Technical Support	31

Required Materials

Introduction

This section identifies the kits, reagents, and equipment needed for this protocol. Vendor and catalog information are provided in the following tables.

SOLiD™ Fragment Library Construction Kit

The following components are needed from the SOLiD[™] Fragment Library Construction Kit (Applied Biosystems 4443473).

End polishing Enzyme 1
End polishing Enzyme 2
5X End-Polishing Buffer
5X T4 Ligase Buffer
T4 Ligase
Platinum® PCR Amplification Mix
SOLiD [™] Library Column Purification Kit with B2-S and B2-L buffers

SOLiD™ Fragment Library Oligos Kit

The following components are needed from the SOLiD™ Fragment Library Oligos Kit (Applied Biosystems 4401151).

P2 adaptor, 50 μM	
Library PCR Primer 1	
Library PCR Primer 2	

SOLiD™ Library TaqMan® Quantitation Kit

All components are needed from the SOLiD[™] Library TaqMan[®] Quantitation Kit (Applied Biosystems 4449639)

Invitrogen Reagents

The following Invitrogen reagents are sold separately.

Product	Catalog no.	
UltraPure [™] DNase/RNase-Free Distilled Water	10977-015	
dATP, 100 mM	10216-018	
dTTP, 100 mM	10219-012	
dGTP, 100 mM	10218-014	
E-Gel® EX Gels, 2% agarose	G4010-02	
E-Gel [®] 50 bp DNA Ladder	10488-099	
Optional if performing the lambda DNA spike-in:		
Lambda DNA	25250-010	
Alu I restriction enzyme	45200-029	
PureLink™ PCR Purification Kit	K3100-01	

Applied Biosystems Reagents

The following Applied Biosystems reagents are sold separately.

Product	Catalog no.
Exo– Klenow Fragment (5 U/μL)	AM2006
0.5 M EDTA, pH 8.0	AM9260G
Hi-Di™ Formamide	4440753
Cells-to-CpG [™] Bisulfite Conversion Kit (50 rxns)	4445249
8-strip PCR Tubes and Caps, RNase-free	AM12230
Low TE buffer	4389764

Additional Reagents

The following reagents from other vendors are sold separately.

Product	Vendor	Catalog no.
5-methyl-dCTP, 100 mM	TriLink (USA)	N-2026
Alternate source of 5-methyl-dCTP	Jena Bioscience (Germany)	NU-1125
Amicon® Ultra-0.5 10K devices	Millipore	UFC501024
Agencourt® AMPure® XP beads	Beckman Coulter	A63880
DNA LoBind Tubes, 1.5 mL	Eppendorf	022431021
Covaris [™] microTUBEs [™] (snap-cap)	Covaris	520045
Isopropanol, 100%	Various	Various
Ethanol, 100%	Various	Various
DTT, molecular biology grade	Various	Various

P1 Adaptor Oligos

The P1 adaptor oligos are two single-stranded oligos that are annealed together to make the double-stranded methyl-P1 adaptor. One of the adaptor oligos (P1A) is methylated; the other (P1B) is not.

The oligos can be manufactured by any supplier capable of providing custom oligos containing 5-methyl-cytidine residues at \geq 90% purity (typically HPLC purified). IDT (Coralville, Iowa, USA) is one such supplier.

The sequences that should be ordered are:

P1A Adaptor Oligo (top strand, 41 bp, 5-methyl-C denoted as "M" below):

5'- MMAMTAMGMMTMMGMTTTMMTMTATGGGMAGTMGGTGAT -3'

P1B Adaptor Oligo (bottom strand, 43 bp, unmethylated)

5´- ATCACCGACTGCCCATAGAGAGGAAAGCGGAGGCGTAGTGGTT -3´

Equipment

The following equipment is required.

Microcentrifuge (e.g., Eppendorf® 5417R tabletop microcentrifuge)

Thermal Cycler (e.g., Applied Biosystems GeneAmp® PCR System 9700)

NanoDrop® ND-1000 Spectrophotometer

Covaris[™] S2 ultrasonicator

Agilent 2100 Bioanalyzer

Vortexer

TaqMan®-compatible qPCR instrument

Magnetic rack (if using Agencourt® AMPure® XP Beads for size selection)

E-Gel[®] iBase[™] Power System

 $E\text{-}Gel^{\mathbb{B}}\ Safe\ Imager^{^{\text{\tiny{TM}}}}\ Real\text{-}Time\ Transilluminator$

Heat block

Description of the Protocol

Protocol Overview

This protocol provides detailed steps to prepare a bisulfite-converted fragment DNA library for sequencing using the SOLiD $^{\text{\tiny{IM}}}$ 4 System. Starting with 5 μ g of purified human genomic DNA, the protocol can yield a library of sufficient complexity for genome-wide measurement of the cytosine methylation pattern at single-nucleotide resolution. This protocol uses products from Applied Biosystems, Invitrogen, and other vendors in a streamlined and validated workflow.

Please refer to the SOLiD[™] 4 System Library Preparation Guide, March 2010, pages 16–40 and associated Appendices for additional information on fragment library preparation.

Workflow Summary

Shear the DNA

To begin, sonicate the input DNA into small fragments with a mean fragment size of 165 bp (range of 150–180 bp) using a Covaris $^{\text{\tiny M}}$ S2 System. Conditions have been tested for shearing 5 μ g DNA in a total volume of 100 μ L. For certain DNA samples, optimizing the shearing protocol may be necessary.

End-Repair the DNA

Next, use End Polishing Enzyme 1 and End Polishing Enzyme 2 to convert DNA that has damaged or incompatible 5'-protruding and/or 3'-protruding ends to 5'-phosphorylated, blunt-ended DNA. For bisulfite-compatible library construction, a dNTP mix lacking dCTP must be used. You can create such a mix by mixing dATP, dTTP, and dGTP each at a final concentration of 10 mM.

Purify the DNA with the SOLiD™ Library Column Purification Kit
PureLink™ columns are supplied with B2-S and B2-L buffers in the SOLiD™
Library Column Purification Kit. This kit is uniquely designed for SOLiD™
workflows and is used at different steps throughout this protocol.

Ligate Methyl-P1 and P2 Adaptors to the DNA

Next, ligate double-stranded P1 and P2 Adaptors to the ends of the end-repaired DNA. The non-methylated P2 adaptor is included in the SOLiD Fragment Library Oligos Kit. For bisulfite-compatible library construction, a modified P1 adaptor containing methylated cytosines must be annealed from high-purity single-stranded oligos purchased separately. After ligation, purify samples using the SOLiD Library Column Purification Kit.

Nick-Translate the DNA

After ligation, the library molecules undergo nick translation with a nucleotide mix that contains 5-methyl-dCTP in place of normal dCTP. This step yields bisulfite-compatible amplifiable library molecules. After nick translation, purify samples with the $SOLiD^{TM}$ Library Column Purification Kit.

Bisulfite Conversion

The Cells-to- $CpG^{\text{\tiny{TM}}}$ Bisulfite Conversion Kit is the preferred set of reagents to use for this step. Add Hi-Di^M Formamide to denature the DNA sample. The temperature and length of time for bisulfite conversion have been optimized to maximize the conversion efficiency and preserve library molecules.

Desulfonation and Purification

Desulfonate the library at room temperature within a centrifugal filtration device. Immediately afterward, the converted DNA is concentrated and exchanged into Low TE buffer within the same device.

Amplify the Library

Amplify the library molecules using at least four cycles of PCR. The bisulfite conversion process generates many uracil residues within the template DNA strands and also leaves the DNA largely single-stranded. Amplification replaces the uracil bases with thymidine bases and returns the library to a double-stranded state.

Size-Select the DNA

To preserve maximum sequence complexity, size-select the ligated, purified DNA to a broad distribution range (100–350 bp) with Agencourt® AMPure® XP beads. After bead-based sizing, the eluate can be directly quantified with the $SOLiD^{TM}$ Library TaqMan® Quantitation Kit.

Quantitate the Library Using qPCR

Quantitate the library using the SOLiD[™] Library TaqMan[®] Quantitation Kit before proceeding to emulsion PCR.

Methods

Before Starting

Tips and Notes

- Start with 5 µg of DNA. **Note:** The amount of DNA loaded onto a PureLink[™] column should be ≤5 µg to ensure maximum yield. If you want to experiment with larger amounts of starting material, be sure to split the sample over multiple PureLink[™] columns during the purification steps.
- Adjust microcentrifuge speeds and times according to the g-forces specified in the protocols. Applied Biosystems recommends the Eppendorf® 5417R tabletop microcentrifuge.
- Prior to loading, dry spin all PureLinkTM columns at $10,000 \times g$ (13,000 rpm) for 1 minute.
- Perform all steps requiring 0.5-mL and 1.5-mL tubes with Eppendorf® LoBind Tubes.
- Thaw reagents on ice before use.

Step 1. Shear the DNA

Introduction

In this step, you shear the DNA using the Covaris^{$^{\text{TM}}$} S2. Sonication conditions have been optimized for shearing the DNA into a mean fragment size of 165 bp (range of 150–180 bp).

Materials Needed

- Covaris[™] S2 ultrasonicator
- Covaris[™] microTUBEs[™]
- LoBind tubes
- Low TE Buffer



- Set the chiller temperature on the Covaris[™] S2 to 2–5°C to ensure that the temperature of the water bath is ~5°C. **Higher temperatures during fragmentation can be detrimental.** The circulated water chiller should be supplemented with 20% ethylene glycol.
- To load and unload the Covaris[™] microTUBE[™] correctly from the microTUBE holder, see the Covaris[™] S2 user manual.

Shearing the DNA

1. Dilute the DNA to $100 \,\mu L$ in 1X Low TE Buffer in a LoBind tube, as follows:

Table 1: DNA Dilution	
Component	Amount
DNA	5 µg
1X Low TE Buffer	Variable
Total Volume	100 μL

- 2. Place a Covaris[™] microTUBE[™] into the loading station on the Covaris[™] S2. Keep the cap on the tube and use a tapered pipette tip to slowly transfer the 100 µL of DNA sample through the pre-split septa. Be careful not to introduce a bubble into the bottom of the tube.
- 3. Fill the Covaris[™] S2 tank with fresh deionized water to level 12 on the graduated fill line label. The water should cover the visible glass part of the tube.
- 4. Shear the DNA using the following Covaris[™] S2 System conditions:
 - Number of Cycles: 6
 - Bath Temperature: 5°C
 - Bath Temperature Limit: 10°C
 - Mode: Frequency sweeping
 - Water Quality Testing Function: Off
 - Duty cycle: **10**%
 - Intensity: 5
 - Cycles/burst: 100
 - Time: 60 seconds
- 5. Remove the Covaris™ microTUBE™ from the loading station. While keeping the snap-cap on, insert a pipette tip through the pre-split septa, then slowly remove the sheared DNA. Transfer the sheared DNA into a new 1.5-mL LoBind tube.

Optional: Spike-In Lambda DNA Digested with Alu I

Introduction

In this optional step, you add endonuclease-digested phage lambda DNA to your sample as a control.

Materials Needed

- Whole lambda DNA
- *Alu* I restriction enzyme
- PureLink[™] PCR Purification Kit
- Low TE Buffer

Purpose of the Lambda DNA

The addition of endonuclease-digested phage lambda DNA is useful for monitoring the conversion efficiency of the bisulfite reaction. Lambda grown in dcm+ bacteria contains only a very small amount of cytidine methylation (on the internal cytosines of CCWGG motifs). Consequently, the frequency with which all other cytosines in lambda-derived sequences are converted to thymidine after treatment with bisulfite and subsequent amplification provides a good measure of the overall conversion efficiency within a complex sample.

Digestion with an enzyme such as *Alu* I is convenient because it produces a large number of fragments with varying sequences that fall within a useful size for such monitoring. Furthermore, since *Alu* I is a blunt-cutter that cuts between G and C of the 5′...AGCT...3′ palindromic motif, the sequences coming from the lambda "spike" can be identified by their ability to be mapped to the lambda genome and by a characteristic CT (converts to "TT") 5′ dinucleotide or AG 3′ dinucleotide.

Creating and Adding the Lambda DNA Standard

To create a lambda DNA standard:

- 1. Digest 50 μg of whole lambda DNA to completion with 100 units of *Alu* I enzyme in a 250 μl volume in an appropriate buffer at 37°C overnight.
- 2. Purify the digested DNA fragments with the PureLink™ PCR Purification Kit
- 3. Measure the concentration of the purified DNA by spectrophotometry.
- 4. Dilute to a working stock concentration of 10 ng/μL with Low TE Buffer.
- 5. For every 1 μ g of starting material, add 10 ng (1 μ L) of *Alu* I-digested lambda DNA directly to the sheared genomic DNA from the previous page (e.g., add 50 ng of digested lambda DNA to 5 μ g of sheared sample DNA).

Step 2. Prepare the Special dNTP Mixes

Introduction

In this step, you prepare the two special dNTP mixes required to construct bisulfite-compatible $SOLiD^{TM}$ fragment libraries.

Materials Needed

- dATP, 100 mM
- dTTP, 100 mM
- dGTP, 100 mM
- 5-methyl-dCTP, 100 mM (TriLink)*
- 10 mM Tris-HCl (pH 7.5), nuclease-free
- Nuclease-free water



*Jena Bioscience (Germany) also supplies 5-methyl-dCTP but as a 10 mM stock solution (catalog no. NU-1125). If you want to use this more dilute stock, use 50 μL per 100 μL final volume and reduce the amount of nuclease-free water from 74 μL to 29 μL in the following protocol.

dNTP Mix without dCTP

Prepare a 100-µL volume of a dNTP mix without dCTP as follows.

Store in sub-aliquots (8 µL needed per library) at -20°C.

Table 2: dNTP Mix without dCTP		
Component	Volume (µL)	
dATP, 100 mM stock	10	
dTTP, 100 mM stock	10	
dGTP, 100 mM stock	10	
10 mM Tris-HCl (pH 7.5), nuclease-free	6	
Nuclease-free water	64	
Total Volume	100	

dNTP Mix with Methyl-dCTP

Prepare a 100- μ L volume of a dNTP mix with 5-methyl-dCTP in place of normal dCTP, as follows.

Store in sub-aliquots (5 μ L needed per library) at –20 $^{\circ}$ C.

Table 3: dNTP Mix with Methyl dCTP		
Component	Volume (µL)	
dATP, 100 mM stock	5	
dTTP, 100 mM stock	5	
dGTP, 100 mM stock	5	
5-methyl-2'-deoxycytidine -5'-triphosphate (5-methyl-dCTP)	5	
100 mM stock (TriLink)		
10 mM Tris-HCl (pH 7.5), nuclease-free	6	
Nuclease-free water	74	
Total Volume	100	

Step 3. Prepare the Methyl-P1 Adaptor

Introduction

In this step, you prepare a modified P1 adaptor that contains 5-methyl cytidine residues in place of normal non-methylated cytidine residues.

Materials Needed

- P1A adaptor oligo (methylated)
- P1B adaptor oligo (unmethylated)
- 5X T4 DNA Ligase Buffer
- Low TE Buffer



5X T4 Ligase Buffer is used as an annealing buffer in the following procedure.

P1 Adaptor Oligos

The P1 adaptor oligos are two complementary single-stranded oligos that are annealed together to make the double-stranded methyl-P1 adaptor. The P1A adaptor oligo is methylated; P1B is not.

The oligos can be manufactured by any supplier capable of providing custom oligos containing 5-methyl-cytidine residues at ≥ 90% purity (typically HPLC purified). IDT (Coralville, Iowa, USA) is one such supplier.

The sequences that should be ordered are:

P1A Adaptor Oligo (top strand, 41 bp, 5-methyl-C denoted as "M" below):

5'- MMAMTAMGMMTMMGMTTTMMTMTATGGGMAGTMGGTGAT -3'

P1B Adaptor Oligo (bottom strand, 43 bp, unmethylated)

5'- ATCACCGACTGCCCATAGAGAGGGAAAGCGGAGGCGTAGTGGTT -3'

Annealing the Adaptor Oligos

Anneal the adaptor oligos as follows:

- 1. Resuspend each P1 adaptor oligo to ~250 μM in Low TE Buffer.
- 2. The extinction coefficient of the methylated strand is $361,400 \, \text{L/(mole-cm)}$ and that of the non-methylated strand is $430,500 \, \text{L/(mole-cm)}$. Measure the exact concentration of the resuspended oligos by UV spectrophotometry
- 3. Dilute each oligo to a final concentration of 125 μ M, and prepare a mixture of the oligo strands as follows.

Table 4: Methyl-P1 Adaptor Annealing Mix	
Component	Volume (µL)
P1A adaptor oligo, 125 μM	40
P1B adaptor oligo, 125 μM	40
5X T4 Ligase Buffer	20
Total Volume	100



After annealing, the proportions above will yield a solution that contains double-stranded methyl-P1 adaptor at a concentration of $50~\mu M$.

Protocol continued on the next page

Protocol continued from the previous page

4. In a thermal cycler such as the Applied Biosystems GeneAmp® PCR System 9700, heat the solution to 95°C then controllably cool it as shown below to permit the complementary oligos to efficiently anneal into double-stranded adaptor molecules.

Table 5: Cooling steps for annealing methyl-	
P1 adaptor in a thermal cycler	
Temperature (°C)	Time (minutes)
95	5
72	5
60	5
50	3
40	3
30	3
20	3
10	3
4	forever

5. After annealing, the double-stranded methyl-P1 adaptor may be handled and stored alongside the P2 adaptor from the SOLiD™ Fragment Library Oligos Kit. As always, be careful to avoid denaturing the annealed adaptors; do not bring them above 37°C.

Step 4. End-Repair the DNA

Introduction

In this step, you end-repair the sheared DNA.

How End-Repair Works

The 5'-to-3' polymerase and the 3'-to-5' exonuclease activities of End Polishing Enzyme 2 generate blunt-ended DNA fragments. End Polishing Enzyme 1 and ATP phosphorylate the 5'-ends of the blunt-ended DNA to enable subsequent ligation.

Materials Needed

- 5X End-Polishing Buffer
- End-Polishing Enzyme 1, 10 U/μL
- End-Polishing Enzyme 2, 5 U/μL
- 1.5-mL LoBind tubes
- Nuclease-free water

End-Repair Mixture

Combine and mix the following components in a 1.5-mL LoBind tube, and incubate at room temperature for 30 minutes.

Table 6: End-Repair Mix	
Component	Volume (µL)
Sheared DNA (5 µg)	100
5X End-Polishing Buffer	40
dNTP mix without dCTP (from page 12)	8
Nuclease-free Water	32
End-Polishing Enzyme 1, 10 U/μL	4
End-Polishing Enzyme 2, 5 U/μL	16
Total Volume	200

Step 5. Purify the End-Repaired DNA

Introduction

In this step, you purify the end-repaired DNA using components in the $SOLiD^{TM}$ Library Column Purification Kit.

Materials Needed

SOLiD[™] Library Column Purification Kit (included with the SOLiD[™] Fragment Library Construction Kit)



The amount of DNA loaded onto a PureLinkTM column should be $\leq 5~\mu g$ to ensure maximum yield. If you are experimenting with larger amounts of starting material, split the sample over multiple PureLinkTM columns during the purification steps.

Purification Procedure

- 1. Before loading the PureLink[™] column, dry spin at 10,000 × g (13,000 rpm) for 1 minute.
- 2. Add 800 μ L (four volumes) of **Binding Buffer (B2-S)** containing 55% isopropanol to the 200 μ L of end-repaired DNA from the previous page. Mix thoroughly.
- 3. Apply ~700 µL of the DNA/binding buffer mixture to the column.
- 4. Let the column stand for 2 minutes at room temperature.
- 5. Centrifuge the column at ≥10,000 × g (13,000 rpm) for 1 minute and discard the flow-through.
- 6. Repeat steps 3–5 to load the rest of the sample onto the column. After discarding the flow-through, place the column back into the same collection tube.
- 7. Add 650 µL of Wash Buffer (W1) to the column.
- 8. Centrifuge the column at ≥10,000 × g (13,000 rpm) for 2 minutes, then discard the flow-through. Repeat to remove any residual wash buffer.
- 9. Air-dry the column for 2 minutes to evaporate any residual alcohol. Insert the column into a clean 1.5-mL LoBind tube.
- 10. Add 50 μ L of Elution Buffer (E1) to the column to elute the DNA, then let the column stand for 2 minutes.
- 11. Centrifuge the column at ≥10,000 × g (13,000 rpm) for 1 minute. **DO NOT DISCARD!** The flow-through contains your sample.
- 12. Add the eluate back onto the column, then let the column stand for 2 minutes. Centrifuge the column once more at $\geq 10,000 \times g$ (13,000 rpm) for 1 minute and collect the flow-through.
- 13. Quantitate the purified DNA by using 2 μ L of the sample on the NanoDrop® ND-1000 Spectrophotometer.



STOPPING POINT! Store the purified DNA in Elution Buffer (E1) at 4°C or proceed directly to **Ligate Methyl-P1 and P2 Adaptors to the DNA**.

Step 6. Ligate Methyl-P1 and P2 Adaptors to the DNA

Introduction

In this step, you ligate the methyl-P1 adaptor and non-methylated P2 adaptor to the DNA.

Materials Needed

- 5X T4 Ligase Buffer
- T4 Ligase
- Methyl-P1 adaptor (double-stranded, from page 13)
- P2 Adaptor (double-stranded, included in the SOLiD™ Fragment Library Oligos Kit)
- 1.5-mL LoBind tube
- Nuclease-free water

Calculating the Amount of Each Adaptor Needed

Before proceeding, calculate the amount of each double-stranded adaptor needed based on the amount of the DNA from the previous purification procedure, using the formulas below.



If DNA fragments were sheared as described on page 9, the average insert size should be approximately 165 bp (before adaptor ligation).

Formulas:

X pmol/μg DNA = 1 μg DNA × (10⁶ pg/ 1 μg) × (1 pmol/660 pg) × (1/average insert size)

Y μL of adaptor needed = # of μg DNA × (X pmol/1 μg DNA) × 15 × (1 μL adaptor needed / 50 pmol)

Example:

For 1 µg of purified end-repaired DNA with an average insert size of 165 bp

X pmol/μg DNA = 1 μg DNA × (10⁶ pg/ 1 μg) × (1 pmol/660 pg) × (1/165) = 9.2 pmol/μg DNA

Y μL adaptor needed = 1 μg DNA × (9.2 pmol/1 μg DNA) × 15 × (1 μL adaptor needed/50 pmol) = **2.76** μL adaptor needed

Ligating the Adaptors

Combine and mix the following components in a 1.5-mL LoBind tube, and incubate at room temperature for 15 minutes.

Table 7: Ligation Mix	
Component	Volume (μL)
Methyl-P1 Adaptor (ds), 50 pmol/μL	Y
P2 Adaptor (ds), 50 pmol/μL	Y
5X T4 Ligase Buffer	40
DNA	48–50
Nuclease-free Water	Variable
T4 Ligase, 5 U/μL	10
Total	200

Step 7. Purify the Adaptor-Ligated DNA

Introduction

In this step, you purify the adaptor-ligated DNA using components in the $SOLiD^{TM}$ Library Column Purification Kit.

Materials Needed

SOLiD[™] Library Column Purification Kit (included with the SOLiD[™] Fragment Library Construction Kit)

Purification Procedure

- 1. Before loading the PureLink[™] column, dry spin at 10,000 × g (13,000 rpm) for 1 minute.
- 2. Add 800 μ L (four volumes) of **Binding Buffer (B2-L)** containing 40% isopropanol to the 200 μ L of adaptor-ligated DNA from the previous page. Mix thoroughly.
- 3. Apply \sim 700 μ L of the DNA/binding buffer mixture to the column.
- 4. Let the column stand for 2 minutes at room temperature.
- 5. Centrifuge the column at ≥10,000 × g (13,000 rpm) for 1 minute and discard the flow-through.
- 6. Repeat steps 3–5 to load the rest of the sample onto the column. After discarding the flow-through, place the column back into the same collection tube.
- 7. Add 650 µL of Wash Buffer (W1) to the column.
- 8. Centrifuge the column at ≥10,000 × g (13,000 rpm) for 2 minutes, then discard the flow-through. Repeat to remove any residual wash buffer.
- 9. Air-dry the column for 2 minutes to evaporate any residual alcohol. Insert the column into a clean 1.5-mL LoBind tube.
- 10. Add 50 μ L of Elution Buffer (E1) to the column to elute the DNA, then let the column stand for 2 minutes.
- 11. Centrifuge the column at ≥10,000 × g (13,000 rpm) for 1 minute. **DO NOT DISCARD THE ELUATE! The flow-through contains your sample.**
- 12. Add the eluate back onto the column, then let the column stand for 2 minutes. Centrifuge the column once more at $\geq 10,000 \times g$ (13,000 rpm) for 1 minute and collect the flow-through.



STOPPING POINT! Store the purified DNA in Elution Buffer (E1) at 4°C or proceed directly to **Nick-Translate the DNA**.

Step 8. Nick-Translate the DNA

Introduction

In this step, you nick-translate the DNA

Materials Needed

- Exo– Klenow Fragment, 5 U/μL (Applied Biosystems)
- 10 mM DTT
- Tris-HCl (pH 8.0)
- MgCl₂
- NaCl
- 0.5 M EDTA, pH 8.0
- 1.5-mL LoBind tube
- Nuclease-free water

Prepare 10X Nick Translation Buffer

Prepare a stock solution of 10X Nick-Translation Buffer (10 μ L needed per reaction). Store at –20°C.

500~mM Tris-HCl (pH 8.0) 100~mM MgCl $_2$ 500~mM NaCl

Nick Translation Procedure

1. Combine and mix the following components in a 1.5-mL LoBind tube:

Table 8: Nick Translation Mix	
Component	Volume (µL)
DNA	48-50
10X Nick Translation Buffer (prepared as above)	10
10 mM DTT	1
dNTP mix with methyl-dCTP (from page 12)	5
Nuclease-free Water	Variable
Exo– Klenow Fragment, 5 U/μL	4
Total	100

- 2. Incubate at 16°C for 1 hour.
- 3. Stop the reaction by adding 4 μ L or 0.5 M EDTA, pH 8.0.

Step 9. Purify the Nick-Translated DNA

Introduction

In this step, you purify the nick-translated DNA using components in the $SOLiD^{TM}$ Library Column Purification Kit.

Materials Needed

SOLiD[™] Library Column Purification Kit (included with the SOLiD[™] Fragment Library Construction Kit)

Purification Procedure

- 1. Before loading the PureLink[™] column, dry spin at 10,000 × g (13,000 rpm) for 1 minute.
- 2. Add 416 μ L (four volumes) of **Binding Buffer (B2-L)** containing 40% isopropanol to the 104 μ L of nick-translated DNA from the previous page. Mix thoroughly.
- 3. Apply the DNA/binding buffer mixture to a PureLink[™] column.
- 4. Let the column stand for 2 minutes at room temperature.
- 5. Centrifuge the column at ≥10,000 × g (13,000 rpm) for 1 minute and discard the flow-through. Place the column back into the same collection tube.
- 6. Add 650 µL of Wash Buffer (W1) to the column.
- 7. Centrifuge the column at ≥10,000 × g (13,000 rpm) for 2 minutes, then discard the flow-through. Repeat to remove any residual wash buffer.
- 8. Air-dry the column for 2 minutes to evaporate any residual alcohol. Insert the column into a clean 1.5-mL LoBind tube.
- 9. Add 50 μL of Elution Buffer (E1) to the column to elute the DNA, then let the column stand for 2 minutes.
- 10. Centrifuge the column at ≥10,000 × g (13,000 rpm) for 1 minute. **DO NOT DISCARD THE ELUATE! The eluate contains your sample.**
- 11. Add the eluate back onto the column, then let the column stand for 2 minutes. Centrifuge the column once more at $\geq 10,000 \times g$ (13,000 rpm) for 1 minute and collect the flow-through.
- 12. Quantitate the purified DNA by using 2 μ L of the sample on the NanoDrop® ND-1000 Spectrophotometer.



STOPPING POINT! Store the purified DNA in Elution Buffer (E1) at 4°C or proceed directly to **Bisulfite Convert the DNA**.

Step 10. Bisulfite Convert the DNA

Introduction

In this step, you use sodium bisulfite to chemically convert the cytosine bases in the DNA to uracil bases in an overnight reaction. Cytosines methylated at the 5 position are resistant to this chemical process, and remain as methylated Cs. Denaturation is promoted by adding Hi-Di^{M} Formamide to the DNA sample.

Materials Needed

- Cells-to-CpG[™] Bisulfite Conversion Kit
- Hi-Di[™] Formamide
- 1.5-mL LoBind tube
- Nuclease-free water



One Cells-to-CpG[™] Bisulfite Conversion Kit provides enough conversion reagent (5 tubes) to convert 5 libraries.

Preparing the Cells-to-CpG™ Conversion Reagent

Prepare complete Cells-to-CpG[™] Conversion Reagent immediately before use:

- 1. In a sterile 1.5-mL tube, mix 26 μ L of Denaturation Reagent with 800 μ L of nuclease-free water.
- 2. Add this mixture to one opaque brown tube of powdered Conversion Reagent and mix well by vortexing. Perform all subsequent steps in this opaque brown tube.
- 3. Add $50 \,\mu\text{L}$ of Conversion Buffer to the tube of resuspended Conversion Reagent, and mix again by vortexing.
- 4. Place the complete Conversion Reagent tube in a 60°C water bath for 30 minutes. Remove and mix the tube every 10 minutes during incubation.
- 5. Following incubation, mix by vortexing 2–3 more times.

Trace amounts of undissolved material may remain; this is normal. Keep the tube protected from light at room temperature until use.



The following reaction requires an 8-hour (e.g., overnight) incubation, and can be held in the thermal cycler at 4°C for several hours afterwards.

Bisulfite Conversion Protocol

- 1. To each purified, nick-translated 50-μL DNA sample from page 20, add 50 μL of Hi-Di™ Formamide.
- 2. Mix thoroughly and heat the mixture at 95°C for 5 minutes to denature the DNA.
- 3. Immediately place the DNA/formamide sample tube on ice for at least 2 minutes.
- 4. After cooling, split the 100 μL DNA/formamide sample into equal 25 μL subsamples in 4 PCR tubes.
- 5. Transfer the complete Conversion Reagent from its opaque brown tube (previous page) to a non-opaque 1.5-mL microcentrifuge tube. Pulse-spin the tube to force any undissolved material to the bottom of the tube.

Note: In the next step, be careful to pipet Conversion Reagent from the top of the tube to avoid transferring any undissolved solids from the bottom.

- 6. Add 125 μ L of complete Conversion Reagent to each 25 μ L DNA/formamide sample. Mix by pipetting up and down.
- 7. Place the tubes in a thermal cycler. Program the cycler to heat at 50°C for 8 hours, then hold the reaction at 4°C for up to 20 hours
- 8. Following incubation, proceed to **Desulfonate and Purify the Converted DNA** on the next page.

Step 11. Desulfonate and Purify the Converted DNA

Introduction

In this step, you desulfonate and purify the bisulfite-converted DNA.

Materials Needed

- Cells-to-CpG[™] Bisulfite Conversion Kit
- Amicon® Ultra-0.5 10K ultrafiltration device (Millipore)
- Low TE Buffer
- Nuclease-free water



IMPORTANT! Dilute the stock Desulfonation Reagent provided in the Cells-to- $CpG^{\mathbb{T}}$ kit with **nuclease-free water**, NOT ethanol. Add 10 mL of nuclease-free water per bottle of stock Desulfonation Reagent.

Dilution with water is required to maximize the yield of bisulfite-converted library DNA by ultrafiltration-based purification.

DO NOT use the water-diluted Desulfonation Reagent for any purpose other than the bisulfite conversion reaction in this protocol. Desulfonation in water is NOT compatible with the standard silica column-based purification described in the Cells-to-CpG™ kit manual.



- When adding solutions to centrifugal filtration devices such as the Amicon® Ultra-0.5 10K, avoid touching the membranes as this can lead to device failure.
- After the 8-hour conversion reaction, some discoloration (light yellow or brown) of the solution may be evident. The color goes away during the course of the desulfonation and purification steps, and has no apparent effect on the quality of the final library DNA.

Desulfonation and Purification Protocol

Desulfonate and purify each 150- μ L bisulfite conversion reaction (sub-sample) separately using an Amicon® Ultra-0.5 10K ultrafiltration device. All steps are at room temperature

- 1. Dilute the stock Desulfonation Reagent provided in the Cells-to- CpG^{T} kit with **nuclease-free water** as described in the **Important** note above.
- 2. Add 200 μ L of nuclease-free water to the Amicon® device, then add a 150- μ L bisulfite-converted DNA sub-sample.
- 3. Spin at $14,000 \times g$ for 10 minutes and measure the flow-through volume (should be ~300 μ L).
- 4. Replace the flow-through with an equal volume of nuclease-free water added onto the column.
- 5. Repeat steps 3–4 one more time.
- 6. Spin at $14,000 \times g$ for 10 minutes and measure the flow-through (~300 µL).
- 7. Replace the flow-through with an equal volume of diluted Desulfonation Reagent.
- 8. Mix well and then let stand for 15 minutes.

Protocol continued on the next page

Continued from the previous page

- 9. Spin at $14,000 \times g$ for 10 minutes and measure the flow-through (~300 µL).
- 10. Replace the flow-through with an equal volume of nuclease-free water.
- 11. Spin at $14,000 \times g$ for 10 minutes and measure the flow-through (~300 µL).
- 12. Replace the flow-through with an equal volume of Low TE Buffer.
- 13. Repeat steps 12–13 one more time.
- 14. Spin at $14,000 \times g$ for 15 minutes and measure the flow-through. This time, ~350 μ L should flow through.
- 15. Invert the device into a clean receiving tube and elute at $1000 \times g$ for 1 minute; the volume that elutes should be between 20 and 40 μ L.
- 16. Re-pool the four sub-samples of bisulfite-converted DNA into a 1.5-mL LoBind tube.
- 17. The eluted, re-pooled, bisulfite-converted, single-stranded DNA library is now ready for PCR amplification.

Step 12. Amplify the DNA

Introduction

In this step, you amplify the bisulfite-converted DNA in a PCR reaction.

The bisulfite conversion process generates many uracil residues within the template DNA strands and also leaves the DNA largely single-stranded. Amplification replaces the uracil bases with thymidine bases and returns the library to a double-stranded state.

Materials Needed

- Platinum® PCR Amplification Mix
- Library PCR Primer 1
- Library PCR Primer 2
- Thermal cycler
- E-Gel® EX 2% Gel
- E-Gel[®] iBase[™] Power System
- E-Gel® Safe Imager™ Real-Time Transilluminator

Determining the Optimal Number of PCR Cycles

To avoid overamplification, determine the minimum number of PCR cycles needed to generate a visible smear of DNA on an agarose gel, as described below.

1. Prepare a test PCR reaction using $1/10^{th}$ of the eluate from the previous page (typically 8–16 μ L).

Table 9: Test PCR Reaction	
Component	Volume µL
Platinum® PCR Amplification Mix	38
Library PCR Primer 1	1
Library PCR Primer 2	1
DNA library sample (eluate from previous page)	8–16 μL

2. Run the PCR program shown in Table 10. Cycle 4 times, then pause the reaction at 4° C to remove 5 μ L to run on a gel. Continue cycling, pausing at 4° C every 2 cycles to withdraw another 5- μ L sample, up to ~12 cycles. Place each sample in a separate LoBind tube.

Table 10: Test PCR Cycling Program			
Stage	Step	Temp (°C)	Time
Holding	Denature	95	5 min
Cycling (collect sample at 4 cycles and	Denature	95	15 sec
at 2-cycle intervals thereafter up to ~12	Anneal	62	15 sec
cycles)	Extend	70	3 min
Holding	Extend	70	5 min
Holding	Hold	4	∞

3. Run the amplified samples on an E-Gel® EX 2% Gel with an E-Gel® iBase™ and E-Gel® Safe Imager™ Real-Time blue-light transilluminator. Determine the minimum number of cycles needed to generate a **faintly visible smear of DNA**. Usually this will be between 4 and 10 cycles. Then prepare and run the following full-scale reaction using that number of cycles.

Amplification Reaction

1. Prepare a master mix based on the volume of eluate from page 24. The volumes below allow for one additional reaction for a no-template negative control:

Table 11: PCR Master Mix		
	Eluate > 100 μL	Eluate ≤ 100 μL
Component	Volume (µL)	Volume µL
	F = Volume of eluate/100	
Platinum® PCR Amplification Mix	380 × F	380
Library PCR Primer 1	10 × F	10
Library PCR Primer 2	10 × F	10
Total	$400 \times F$	400

- 2. If the volume of the eluate is:
 - $\leq 100~\mu L$, add $400~\mu L$ of master mix to the eluate, then distribute equally into four PCR reaction tubes.
 - $> 100~\mu L$, add $400~\mu L$ of master mix to every $100~\mu L$ of eluate, then distribute equally in 100– $125~\mu L$ aliquots to multiple PCR reaction tubes.
- 3. Prepare and run the following PCR program, using the number of cycles determined previously.

Table 12: PCR Cycling Program				
Stage	Step	Temp (°C)	Time	
Holding	Denature	95	5 min	
Cycling (number of	Denature	95	15 sec	
cycles determined by	Anneal	62	15 sec	
the previous test PCR)	Extend	70	3 min	
Holding	Extend	70	5 min	
Holding	Hold	4	∞	

4. After cycling, pool the reactions into 200–250 μL volumes in multiple 1.5-mL LoBind tubes.

Step 13. Purify the Amplified DNA

Introduction

In this step, you purify the amplified DNA using components in the SOLiD[™] Library Column Purification Kit.


Materials Needed

SOLiD[™] Library Column Purification Kit (included with the SOLiD[™] Fragment Library Construction Kit)

Purification Procedure

Perform the following workflow for each 200–250 µL pooled PCR volume from the previous page.

- Before loading the PureLink[™] columns, dry spin at 10,000 × g (13,000 rpm) 1. for 1 minute.
- Add four volumes of **Binding Buffer (B2-L)** containing 40% isopropanol to each 200–250 µL volume of amplified DNA from the previous page. Mix thoroughly.
- For each sample, apply ~700 µL of the DNA/binding buffer mixture to a PureLink[™] column.
- 4. Let the columns stand for 2 minutes at room temperature.
- Centrifuge the columns at $\geq 10,000 \times g$ (13,000 rpm) for 1 minute and 5. discard the flow-through.
- Repeat steps 3–5 to load the rest of each sample onto the column. Place the column back into the same collection tube.
- 7. Add 650 µL of Wash Buffer (W1) to each column.
- Centrifuge the columns at $\geq 10,000 \times g$ (13,000 rpm) for 2 minutes, then discard the flow-through. Repeat to remove any residual wash buffer.
- 9. Air-dry the columns for 2 minutes to evaporate any residual alcohol. Insert each column into a clean 1.5-mL LoBind tube.
- 10. Add 50 µL of Elution Buffer (E1) to each column to elute the DNA, then let the column stand for 2 minutes.
- 11. Centrifuge the columns at ≥10,000 × g (13,000 rpm) for 1 minute. **DO NOT** DISCARD THE ELUATE! The eluate contains your sample.
- 12. Add the eluate back onto each column, then let the columns stand for 2 minutes. Centrifuge the columns once more at ≥10,000 × g (13,000 rpm) for 1 minute and collect the flow-through.
- 13. Pool the eluted DNA from the PureLink[™] columns into a LoBind tube.
- 14. Quantitate the purified DNA library by using 2 µL of the sample on the NanoDrop® ND-1000 Spectrophotometer.

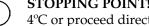

STOPPING POINT! Store the purified DNA in Elution Buffer (E1) at 4°C or proceed directly to **Size-Select the DNA**.

Step 14. Size-Select the DNA

Introduction

In this step, you size-select the DNA.

Materials Needed

- AMPure® XP Beads
- 1.5-mL LoBind tubes
- Magnetic rack
- 70% ethanol, freshly prepared
- Low TE Buffer

Bead-Based Selection Using AMPure® XP Beads

This method selects for amplified adaptor-ligated DNA molecules that are between ~100 and ~350 bp in length, with an average length of~200 bp.

- 1. Add 0.8 volumes (relative to the volume of DNA solution) of Agencourt AMPure® XP beads to the sample. For example, if the pooled, eluted DNA from the previous page is in 100 μ L, add 80 μ L of bead solution.
- 2. Mix well and incubate the solution for 5 minutes at room temperature.
- 3. Place the tube of beads in a magnetic rack to separate the beads from the solution. Wait 2 minutes for the solution to clear before proceeding to the next step.
- 4. Keep the tube in place in the magnetic rack with the beads held against the tube wall and carefully remove and RETAIN the supernatant in a clean 1.5-mL LoBind tube. The supernatant contains your sample.
- 5. To this supernatant, add 1 volume (original volume) of Agencourt AMPure® XP beads (100 μL in the example from step 1 above).
- 6. Mix well and incubate the beads and DNA solution for 5 minutes at room temperature.
- 7. Place the tube of beads in a magnetic rack to separate the beads from the solution. Wait 2 minutes for the solution to clear before proceeding to the next step.
- 8. Keep the tube in place in the magnetic rack with the beads held against the tube wall and gently remove and DISCARD the supernatant.
- 9. With the tube still in place in the magnetic rack, gently add 200 µL of freshly prepared 70% ethanol. Keeping the tube in place on the magnet, let the bead pellet soak in the ethanol solution for 2 minutes at room temperature. Gently aspirate out and discard the ethanol solution.
- 10. Repeat this washing procedure twice with clean 70% ethanol each time.
- 11. Keeping the tube in place on the magnet, air-dry the beads (with the tube cap open) for 5 minutes at room temperature.
- 12. Elute the DNA by adding $22 \,\mu\text{L}$ of Low TE Buffer, vortexing for $10 \,\text{seconds}$, pulse-spinning the solution to the bottom of the tube, and ensuring homogeneity by pipetting the solution up and down several times.

Protocol continued on the next page

Protocol continued from the previous page

- 13. Place the tube of beads in the magnetic rack to separate the beads from the solution. Wait 2 minutes for the solution to clear before proceeding to the next step.
- 14. Carefully transfer the supernatant to a new 1.5-mL LoBind tube.
- 15. Place the eluted sample again in the magnetic rack to separate any remaining beads from the solution. Wait for the solution to clear before proceeding to the next step.
- 16. Carefully transfer the supernatant to a new 1.5-mL LoBind tube. The library is now ready for quantitation.



STOPPING POINT! Store the purified DNA at 4°C or proceed directly to **Quantitate the Library**.

Step 15. Quantitate the Library

SOLiD™ Library TaqMan® Quantitation Kit

For accurate library quantitation, quantitative PCR (qPCR) is strongly recommended. The SOLiD™ Library TaqMan® Quantitation Kit is available from Applied Biosystems. See the user manual for that kit for instructions.



STOPPING POINT! Following quantitation, store the purified DNA at −20°C or proceed directly to emulsion PCR as described in the Applied Biosystems SOLiD[™] 4 System Templated Bead Preparation Guide (PN 4448378) or the Applied Biosystems SOLiD[™] EZ Bead[™] Emulsifier Getting Started Guide (PN 4441486).

Technical Support

Support Information

For the latest services and support information for all locations, go to: www.appliedbiosystems.com

At the Applied Biosystems web site, you can:

- Access worldwide telephone and fax numbers to contact Applied Biosystems Technical Support and Sales facilities.
- Search through frequently asked questions (FAQs).
- Submit a question directly to Technical Support.
- Order Applied Biosystems user documents, SDSs, certificates of analysis, and other related documents.
- Download PDF documents.
- Obtain information about customer training.
- Download software updates and patches.

Part number: MAN0003271 Revision date: 2 December 2010



International Sales