



## HDM97 *Pocket* Reference Meter Series



## User Manual

**HDM97BL**  
**HDM97BM**  
**HDM97BN**  
**HDM97BO**  
**HDM97BP**  
**HDM97BQ**

# Preface

The information contained in this is subject to change without notice. IBP Medical GmbH, its distributors and subsidiaries take no responsibility for any errors or omissions in this document. The contained software is being delivered on the basis of a general licence contract or in single license. Use or reproduction of the software is allowed only in agreement with the contractual arrangements. Whoever transfers this software and/or this manual on magnetic tape, diskette or any other media, except for the purpose of own use, without written authorization of the IBP Medical GmbH, is liable to prosecution.

Copyright (C) 2009 IBP Medical GmbH.  
All rights reserved

All brand names mentioned in this document are property of their respective owners.

Publishers:	IBP Medical GmbH Ikarusallee 15 D 30179 Hannover Germany	IBP Medical Inc 4113 W. St. Charles Ave. Phoenix, AZ 85041 USA
Phone:	+49 511 651647	1-865-686-8650
Fax:	+49 511 652284	1-866-243-0187
Internet:	<a href="http://www.ibpmedical.com">http://www.ibpmedical.com</a>	<a href="http://www.ibpmedical.com">http://www.ibpmedical.com</a>
E-Mail:	<a href="mailto:support@ibpmedical.com">support@ibpmedical.com</a>	<a href="mailto:america@ibpmedical.com">america@ibpmedical.com</a>
Manual authors:	Dipl. Ing. Werner Pfingstmann Holger Kleinert	
Document Number	DHDM97.0220.01	
Revision history	Rev. 0. 29.05.2008 Rev. 1. 18.03.2009 Update Hard- und Software release 2.0 Rev. 2. 01.10.2009 Update Flow sensor Rev. 3. 04.11.2009 Update Flowsensor Cleaning	

**This manual is only valid for the software revisions listed below:**

<b>HDM97BL</b>	<b>V 2.00 ...2.19</b>
<b>HDM97BM</b>	<b>V 2.00 ...2.19</b>
<b>HDM97BN</b>	<b>V 2.00 ...2.19</b>
<b>HDM97BO</b>	<b>V 2.00 ...2.19</b>
<b>HDM97BP</b>	<b>V 2.00 ...2.19</b>
<b>HDM97BO</b>	<b>V 2.00 ...2.19</b>
<b>HDM97BQ</b>	<b>V 2.00 ...2.19</b>

<b>IBPView</b>	<b>V 2.00</b>
----------------	---------------







# SAFETY INSTRUCTIONS



---

For your and your patients safety read and consider the safety instructions below carefully

---

- Read the entire manual carefully before using the *HDM97Pocket* .
- Keep the instrument away from unauthorised persons.
- Only use the *HDM97Pocket* as reference instrument to measure Conductivity, Temperature, Pressure, Flow and pH depending to it's equipped features.
- Never use the *HDM97Pocket* as a replacement of sensors of a medical device.
- Do not use the *HDM97Pocket* in conjunction with a medical device during a treatment.
- Only use the instrument in a dry environment and do not touch it with wet hands.
- Make sure that no liquids intrude inside the instrument or it's inlets / outlets.
- Always use a clean protection filter for the pressure transducer inlet.
- Prevent every mechanical overstraining of electrical wires. Never buckle or pull the cable of the conductivity probe.
- If the acquired values seem to be not believable, make sure that the *HDM97Pocket* is not defective
- Prevent electrostatic discharge on the connectors. This can lead to substantial damage of the instrument. Make sure to be completely discharged before touching the connectors or cables connected to the instrument.
- Adjust the meter only, if you are familiar with the consequences of the adjustment. Consider the hints for reference solutions in the user manual.
- Potentials above 42V against earth ground are dangerous. This potentials can lead to electrical shock and therefore to health hazards. Make sure that none of the connectors has higher voltage than mentioned before.
- The direct connection of earth potential to any connector of the instrument is not applicable for safety purposes and may substantially damage the instrument. Make sure that the medical device itself is properly connected to earth ground according to its instruction manual.
- Make sure that the instrument never heats over 60°C / 140°F. Prevent direct sunlight.
- Never sterilize the instrument using an autoclave. Danger of explosion of the battery!
- Never throw the meter into a fire. Danger of explosion of the battery!
- Never open the meter. There are no parts inside you can repair.
- Never try to replace or repair the internal battery. Handled in the wrong way these part is dangerous. Extreme risk of fire and explosion!

**Table of content**

<b>Introduction .....</b>	<b>2</b>
<b>Application .....</b>	<b>2</b>
<b>EC DECLARATION OF CONFORMITY.....</b>	<b>3</b>
<b>Delivery Content .....</b>	<b>4</b>
<b>Warranty .....</b>	<b>5</b>
<b>Product Overview .....</b>	<b>6</b>
The versions at a glance .....	6
Measuring Channels .....	7
Charging the battery .....	8
Performing measurements while the battery is being charged .....	8
Flow-through adapter for conductivity and temperature measurement .....	8
Keypad .....	9
Switching on the instrument .....	10
Display test during start up.....	10
Power saving function .....	10
Displaying battery charge progress and finish .....	10
Battery gauge display during operation.....	11
Battery aging .....	11
Battery replacement – ONLY BY MANUFACTURER .....	11
Value adjustment.....	14
<b>Measuring .....</b>	<b>15</b>
Conductivity and temperature measurement .....	15
Fundamentals of temperature compensation.....	15
Appliance of temperature coefficients .....	15
Temperature compensation in dialysis.....	17
Temperature measurement.....	19
Pressure measurement .....	19
Pressure drop measurement.....	20
pH measurement.....	20
Fundamental information about pH measurement.....	20
<b>Stopwatch (TIMER) .....</b>	<b>21</b>
Flow measurement.....	21
<b>Attenuation of reading.....</b>	<b>22</b>
<b>Adjustment and Calibration .....</b>	<b>23</b>
<b>Adjustment and Calibration .....</b>	<b>23</b>
Precautions .....	23
Definitions.....	23
General.....	23
Handling of Reference Solutions.....	24
Adjusting the HDM97 Pocket .....	24
<b>Maintenance of HDM97 Pocket.....</b>	<b>27</b>
<b>Troubleshooting.....</b>	<b>30</b>
<b>Technical data .....</b>	<b>31</b>
<b>PC Interface .....</b>	<b>33</b>
USB Interface .....	33
Interface parameters for serial COM port emulation .....	33
Driver installation, example for <i>Microsoft® Windows® XP Service Pack 2</i> .....	34
Example.....	36
<b>IBPView.....</b>	<b>37</b>
What is IBPView .....	37
Installation .....	38
Quickstart .....	38
Disconnect USB .....	38
Main Window .....	38
Selecting hardware.....	39
Assigning graphs to hardware channels .....	40
List of channels .....	41

## Introduction

The HDM97 *Pocket* meter series are reference instrument designed for the calibration of medical devices such as hemodialysis machines. Please read this entire manual carefully before the first use of the instrument.

## Application

Depending on the version the HDM97 *Pocket* may be equipped with up to 5 acquisition channels:

Channel	HDM97BL	HDM97BM	HDM97BN	HDM97BO	HDM97BP	HDM97BQ
Conductivity			✓	✓	✓	✓
Temperature			✓	✓	✓	✓
Timer	✓	✓	✓	✓	✓	✓
Pressure		✓		✓	✓	✓
pH					✓	
Flow	✓					✓

The instrument was designed to be used as a dialysis machine maintenance instrument, but can also be used as laboratory measuring instrument as well as instrument for environmental purposes.

This instrument displays measuring values numeric. Graphical displays are available through a PC or Notebook using the optional Software *IBPView*.

Using the USB communication interface it is possible to obtain measuring values regardless of the instrument display's content. These values can be displayed in any way or stored for further processing.

***EC DECLARATION OF CONFORMITY***  
*according to the Council Directive 93/42/EEC concerning medical devices*

**Manufacturer**

IBP Medical GmbH  
Ikarusallee 15  
30179 Hannover

*declare under our sole responsibility that the following products:*

***Dialysis Reference Meter***

**Version** ***HDM97BL***  
***HDM97BM***  
***HDM97BN***  
***HDM97BO***  
***HDM97BP***  
***HDM97BQ***  
***HDM99XP***

*Product Code: 31.0XXX.XX*

*X = Depending on Version*

**meet the provisions of the Council Directive 93/42/EEC concerning medical devices which apply to them.**

***Product Classification***

II B

**Notified Body**

Medcert GmbH  
Vorsetzen 35  
20459 Hamburg

***Identification Number***

0482

Hannover 18.04.09

Responsible:  
Dipl. Ing. Werner Pfingstmann

## Delivery Content

Pos.	Description	Version	Standard	Option
1	HDM97 <i>Pocket</i> meter <i>BL, BM, BN, BO, BP, BQ</i>		✓	
2	Conductivity/Temperature Probe	<i>BN, BO, BP, BQ</i>	✓	
3	Dialyser Line Adapter	<i>BN, BO, BP, BQ</i>	✓	
4	Hose for pressure measurement	<i>BM, BO, BP, BQ</i>	✓	
5	14,00 mS/cm Conductivity Reference Solution	<i>BN, BO, BP, BQ</i>	✓	
6	Battery charger with wall plug adapters	All	✓	
7	USB interface cable	All	✓	
8	CD - HDM97 <i>Pocket</i> with USB drivers, Software and User Manual	All	✓	
9	Safety instructions and quick start guide	All	✓	
10	Flow Sensor	<i>BL</i>	✓	
11	Flow Sensor	<i>BQ</i>		✓
12	pH Probe with buffer solution pH6, 7 and 8	<i>BP</i>		✓
13	Carrying case	All		✓
14	Holder Infusion pole	All		✓
15	CD - IBPView Data Acquisition Software	All		✓

## Warranty

IBP Medical GmbH warrants that it will repair or replace, at its option, any defective or malfunctioning part without charge for the terms listed below. Parts used for repair or replacement are warranted for the remaining warranty period only.

The user must deliver, at its own expense, the product to IBP Medical GmbH.

<b>Parts</b>	<b>Warranty –Period*</b>
HDM97 <i>Pocket</i>	24 Month
pH-Probe	6 Month

\* from date of delivery note

### The warranty does not cover:

- Disposable parts as battery or pH electrode
- Annually calibration
- Cell cleaning
- Defects caused by:
  1. Modification, alteration, repair or service of the product by anyone other than IBP Medical or an authorized service center
  2. Misuse due to negligence or accident
  3. Operation or maintenance of the product in a manner contrary to IBP instructions

Any express warranty not provided herein, and any remedy for breach of contract that but for this provision might arise by implication or operation of law, is hereby excluded and disclaimed. The implied warranties of merchantability and of fitness for any particular purpose are expressly limited to the terms mentioned above. Some states do not allow limitations on the duration of an implied warranty, so the above limitation may not apply to you.

Under no circumstances shall IBP Medical GmbH be liable to the original purchaser or to any other person for any special or consequential damages, whether arising out of breach of warranty, breach of contract, or otherwise. Some states do not allow the exclusion or limitation of special or consequential damages, so the above exclusion or limitation may not apply to you.

This warranty gives you specific legal rights, and you may also have other rights that vary from state to state.

For further warranty information, contact IBP Medical GmbH.

## Product Overview

### HDM97<sub>BL</sub>



Flow

### HDM97<sub>BM</sub>



Pressure

### HDM97<sub>BN</sub>



Conductivity  
Temperature

### HDM97<sub>BO</sub>



Pressure Conductivity  
Temperature

### HDM97<sub>BP</sub>



pH Pressure Conductivity  
Temperature

### HDM97<sub>BQ</sub>



Flow Pressure Conductivity  
Temperature

The versions at a glance

### HDM97<sub>BL</sub>

This version is suited for usage with a HDU FLOW Sensor and is also equipped with a timer function

### HDM97<sub>BM</sub>

This version is equipped with a high resolution pressure transducer, has a pressure drop measurement function and a timer function. The high precision pressure transducer is not waterproof!

### HDM97<sub>BN</sub>

This version is equipped with the conductivity / temperature probe only. It also has the timer function built in.

### HDM97<sub>BO</sub>

This version is equipped with the conductivity / temperature probe and a standard precision water proof pressure transducer. It also has the pressure drop measurement and the timer function built in.

### HDM97<sub>BP</sub>

This version is equipped with the conductivity / temperature probe and a standard precision water proof pressure transducer. It also has the pressure drop measurement and the timer function built in. It can also measure pH with standard pH sensors.

### HDM97<sub>BQ</sub>

This version is equipped with the conductivity / temperature probe and a standard precision water proof pressure transducer. The high precision pressure transducer is available as "Option H" for this version. This high precision pressure transducer is not waterproof. The meter can also measure flow by using the HDU FLOW probe. It also has the pressure drop measurement and the timer function built in.

## Measuring Channels

Only use accessories provided with the HDM97 *Pocket*. Use of other devices may damage the instrument and will void the warranty.

### ***Conductivity/Temperature Probe***

The conductivity/temperature transducer is a quadropole design for greater accuracy and longevity than other designs.

### ***Standard precision pressure transducer***

This pressure transducer is watertight. However it is recommended to keep fluids away from the sensor because residues like salt can falsify measurement results. In this case the sensor needs to be flushed with deionized water.

### ***High precision pressure transducer***

To avoid damage or inaccurate measurement, ensure that no fluids enter the instrument. We recommend the use of a Transducer Protector for this purpose. Never flush the sensor.

### ***Pressure drop measurement***

This is not a real channel, it is a function for measuring pressure loss in systems for a specified time.

### ***Timer***

This is not a real channel, it is a function to help performing anything where time is a crucial element.

### ***pH***

The input connector is designed exclusively for pH electrodes. It is optimized for glass electrode single-rod measuring cells with a slope of  $\Delta E$  of 59mV at 20°C. Its absolute maximum allowable input voltage span extends from -3 Volts to +3 Volts with respect to BNC Shield. Applying any voltage beyond these boundaries will possibly destroy the instrument.

### ***Flow***

This connection is designed to be used with the HDU FLOW probe only. Do not connect any other devices to this port even when the connector may fit.

## Charging the battery

The HDM97 *Pocket* is internally equipped with a high performance Lithium-Ion-Polymer battery. It can be charged using the external power supply, a PC or a laptop which is connected via the USB cable to the instrument. The battery of the HDM97 *Pocket* can only be charged from USB ports which are capable of delivering 500 mA of current.



This is appropriate to most connectors of laptops or PC's. Connecting the meter to a passive (e.g. not externally powered) USB hub, the battery can not be charged and the communication may not work well. During the charging process the battery symbol in the display is animated.

## Performing measurements while the battery is being charged

The HDM97 *Pocket* can be used for measuring while the battery is being charged. To maintain the dialysis machine and the patient's safety the USB part is isolated to the appliance part of the meter withstanding a voltage of 4000 Volts AC or 6000 Volts DC for at least one minute according to EN60601-1. The appliance part consists of the conductivity, temperature, pressure and pH channel and its connectors, depending on the version.



**ATTENTION: The USB connector is not part of the appliance part.**

## Flow-through adapter for conductivity and temperature measurement



The adapter is suitable for measuring pressure, temperature and conductivity in flow-through mode. With immersion usage of the probe it is suitable for measuring temperature and conductivity concurrently. The preferred orientation of the adapter is vertical. The axial connector must point down (see picture). In flow-through mode the axial connector is to be used as inlet meanwhile the side connector is to be used as outlet. The connection to the pressure transducer inside the HDM97 *Pocket* is accomplished by the lateral connection above the inlet. Before measurement it is necessary to shake the adapter slightly, so that air bubbles trapped inside can pass off through the outlet.

For immersion usage the probe must be unscrewed from the flow-through adapter. To accomplish that do not turn the upper shank, the probe may be damaged. Only turn the nut as shown in the picture left.

**Never touch the electrode surfaces with your fingers or other objects.**

**Do not buckle the cable of the probe. The minimum allowed bending radius is 3 cm e.g. 1¼ inches.**

## Keypad

**Main functions**

The keypads main functions of the HDM97 *Pocket* are listed below:

- Key  for switching the instrument on and off
- Key  for selecting the channel CD/TP ; TP ; PRESSURE ; pH
- Key  for selecting the unit to display.
- Key  leads to channel dependent special functions
- Key  to tare the pressure transducer

The keypad layout is shown in the left picture.

The keys  ,  and  have different sub functions depending on the channel currently displayed. These sub functions are explained in chapter "Measuring"

**Additional functions**

The yellow keys below the display can be used to set values and navigate through menu options or setting. These functions become available when the symbols shown in the right picture appear at the bottom of the display.






Cancels the current setting or selection and returns to measuring mode.



Decreases a value or selects the previous option in a menu or a setting.



Increases a value or selects the next option in a menu or a setting.



Applies a value or a selects an option in a menu or a setting.



## Battery gauge display during operation

If nothing is connected to the USB port during operation the internal battery is used to supply power to the instrument. The battery symbol in the left top corner of the display shows the remaining capacity of the internal battery. The combinations have the meanings listed below:



The battery is 100% full.



The battery has approx. 2/3 of its capacity left.



The battery has approx. 1/3 of its capacity left.



The battery is empty, approx 10% is left.



If this symbol starts flashing the battery is totally empty and the meter will switch off shortly.

## Battery aging

The battery technology of the HDM97 **Pocket** is based on Lithium-Ion-Polymer technology. This technology has many advantages compared with Nickel-Metal-Hydride technology which are lower weight, smaller size, longer life, no memory effect and higher efficiency. However, after a couple of years (5 years estimated) the battery's capacity will start decreasing.

## Battery replacement – ONLY BY MANUFACTURER

Do not try to replace the battery by yourself. The battery is mounted directly onto the main PCB inside the HDM97 **Pocket**, and the main PCB itself contains the battery protection circuit. If the battery is not handled in the right way, there is extremely danger of fire and explosion. Therefore, this battery can only be replaced from IBP personnel who have run through a special training in handling Lithium-Ion batteries.

## Basic settings

To invoke the installation menu please follow the steps below:

1. Switch off the meter using the  key.
2. Switch it on again using the same  key.
3. Immediately press the  key once during the display test.
4. After this procedure the instruments starts the basic settings dialog instead of entering any measurement functions.
5. In the basic settings dialog the user is able to adjust a variety of settings.  
These are:

Parameter	Meter Versions	Description
AUTO-OFF	All	Set the delay time for the auto power off function or disables it.
BEEP	All	enables / disables the keypad acknowledge beep
CAL-RES	All	Overwrites the user calibration dataset with the device initial calibration dataset.
LANGUAGE	All	Adjusts the display language.
PHT	<i>BM,BO,BP,BQ</i>	Target time adjustment for pressure hold test (PHT)
TIMER	All	enables / disables the stopwatch function
Flow	<i>BL, BQ</i>	enables / disables the Flow channel
pH	<i>BP</i>	enables / disables the pH channel
TEMP	<i>BN, BO, BP, BQ</i>	enables / disables the sole readout of the temperature channel

▼ → **AUTO-OFF**

	→		OFF	→	disables function
	→		30 MIN	→	30 minutes delay
	→		60 MIN	→	60 minutes delay
	→		90 MIN	→	90 minutes delay

Then  to apply and save this setting,  
or  to discard this setting.

▼ → **BEEP**

	→		ON	→	keypad beep enabled
	→		OFF	→	keypad beep disabled

Then  to apply and save this setting,  
or  to discard this setting.

- ▼ → **CAL-RES**  
 → Press 5 seconds until the counter in the display reaches zero. Then a calibration dataset reset will be performed. This takes a couple of seconds. The instrument performs a restart afterwards.
- ▼ → **LANGUAGE**  
 → ▲ DEUTSCH → german language  
 → ▼ ENGLISH → english language  
 Then  to apply save this setting,  
 or  to discard this setting.
- ▼ → **TIMER**  
 → ▲ ON → stopwatch function enabled  
 → ▼ OFF → stopwatch function disabled  
 Then  to apply and save this setting,  
 or  to discard this setting.
- ▼ → **PHT (pressure hold test)**  
 → ▲ increases the value  
 → ▼ lowers the value  
 Then  to apply and save this setting,  
 or  to discard this setting.
- ▼ → **PH**  
 → ▲ ON → pH enabled  
 → ▼ OFF → pH disabled  
 Then  to apply save this setting,  
 or  to discard this setting.
- ▼ → **FLOW** only available with version *BQ*  
 → ▲ ON → FLOW enabled  
 → ▼ OFF → FLOW disabled  
 disabled  
 Then  to apply save this setting,  
 or  to discard this setting.
- ▼ → **TEMP**  
 → ▲ ON → sole readout of temperature enabled  
 → ▼ OFF → sole readout of temperature disabled  
 Then  to apply and save this setting,  
 or  to discard this setting.

▼ → **EXIT**

A save dialog will appear, if changes have been made.

→ **SAVE**

→ ▲ YES to apply and save all settings by exiting  
 → ▲ NO to discard all settings by exiting

Then  to exit the basic settings dialog,  
 or  to return to the basic settings dialog.

## Value adjustment

The adjustment of values is done via the navigation keys ▲ and ▼. By pressing and holding down one of these keys the last (right) digit of the value will be incremented or decremented depending on the key. If this digit reaches the boundary it remains 0 and the next digit will begin to change. If the key is being released and pressed again the procedure will start again beginning with the last digit.

### Example showing a value, which has an allowable range from 0.00 to 7.50:

Display: 0.00 ; navigationkey ▲ is hold down:

Order of value incements (every change follows with a delay of 0.3 seconds):

0.01→0.02→0.03→0.04→0.05→0.06→0.07→0.08→0.09 (carry-over)  
 0.10→0.20→0.30→0.40→0.50→0.60→0.70→0.80→0.90 (carry-over)  
 1.00→2.00→3.00→4.00→5.00→6.00→7.00→7.50 (Max value reached)

If the value will exceed the range limits (in this case from 7.00 to 8.00) the value will be truncated to the maximum allowed value, 7.50 in this particular example. The Symbol ▲ will become blank and the meter does not react on the corresponding navigation key. The values can only be changed down to 0.00 using the other ▼ key.

## Measuring

By pressing the key  the user can select between all available channels of the HDM97 *Pocket*. These channels are

- |                                |                   |   |
|--------------------------------|-------------------|---|
| • Conductivity and temperature | Display "CD/TEMP" | only Version <i>BN, BO, BP, BQ</i>            |
| • Temperature                  | Display "TEMP"    | only Version <i>BN, BO, BP, BQ</i> if enabled |
| • Pressure                     | Display "PRES"    | only Version <i>DA, BO, BP, BQ</i>            |
| • pH                           | Display "PH"      | only HDM97 <i>BP</i> , if enabled             |
| • Flow                         | Display "FLOW"    | only HDM97 <i>BQ</i> , if enabled             |
| • Timer                        | Display "Time"    | if enabled                                    |

### Conductivity and temperature measurement

In this mode the display shows "CD/TEMP" in the upper line, the temperature value in the middle line and the conductivity value in the bottom line.

By pressing the key  the user can choose the temperature coefficient.

### Fundamentals of temperature compensation

#### Definition

#### Electrical Conductivity

Electrical conductivity or specific conductivity is a measure of a solution's ability to conduct an electric current. Electrical conductivity of electrolytic solutions is strongly dependent on temperature. In order to compare electrical conductivity measurements at different temperatures, they are standardized to a common reference temperature. This common temperature is 25°C. The dimension unit commonly used is "mS/cm at 25°C"

#### Common temperature

The common temperature is the temperature at which non compensated conductivity measures and compensated conductivity measures have the same value.

#### Temperature coefficient

The temperature coefficient is the relative change of the solution's conductivity depending on it's temperature standardized to a common reference temperature. The dimension unit of this value is measured in %/K.

### Appliance of temperature coefficients

#### **This is a topic that is frequently misunderstood and often neglected in hemodialysis.**

A solution's conductivity will change according to temperature. With increasing temperatures, the measured solution's conductivity will increase, too. To achieve meaningful measurement results, the conductivity value displayed is compensated to 25°C. In other words, the display is always converted to a solution temperature of 25°C. The temperature coefficient which the displayed value is compensated with is expressed as %/°C. Unfortunately however, different solutions also have different temperature coefficients. To achieve an exact display, the instrument will have to be adjusted to the temperature coefficient of the current solution. The commonly used temperature coefficient for dialysates is 2.07 %/K. For naturally occurring solutions, a value of 1.97 %/°C is frequently used. Many measuring devices not specially tailored to dialysis will use this value.

## Temperature compensation

### Temperature compensation using a single temperature coefficient

A single temperature coefficient is used for calculating the conductivity value at the common reference temperature from the actual measured conductivity and the actual measured temperature.

**In order to use one single temperature coefficient the values for 25°C and 37°C must be the same.**

### Dynamic temperature compensation

Dynamic temperature compensation is achieved by calculating a temporary temperature coefficient for the measured temperature using a set of two coefficients for 25°C and 37°C respective. This temporary temperature coefficient is then applied to the actual measured conductivity to calculate the conductivity at the common reference temperature. The following table shows the usage of dynamic temperature coefficients applied to a NaCl solution of 14mS/cm.

Temperature	Conductivity NaCl 14 mS/cm without compensation	Temperature coefficient	Conductivity Readout
25 °C	14,00 mS/cm	<b>2,000 %/K</b>	14,00 mS/cm
31 °C	15,71 mS/cm	2,035 %/K	14,00 mS/cm
37 °C	17,48 mS/cm	<b>2,070 %/K</b>	14,00 mS/cm

**The values 2.00 %/K at 25°C and 2.07 %/K at 37°C are defaults for the respective dialysis machine or values entered by the user. For demonstration purposes the temperature coefficient for 31°C is mentioned, which is calculated by the HDM97Pocket from the values for 25°C and 37°C by linear interpolation.**

## Temperature compensation in dialysis

Listed below are the temperature coefficients used by most major hemodialysis machine manufacturers. We recommend that you double check this data with documentation from your machine's manufacturer

Display	Maschine Manufacturer	Temperature coefficient at 25°C	Temperature coefficient at 37°C
BAXTER	Baxter (Europe Machine)	2,200 %/K	2,200 %/K
BELLCO	Bellco	2,000 %/K	2,100 %/K
B.BRAUN	B.Braun	2,000 %/K	2,100 %/K
FRESENI	Fresenius	2,000 %/K	2,100 %/K
GAMBRO	Gambro (AK-Type Machines)	2,000 %/K	2,070 %/K
GAMBRO	Gambro Phoenix	2,000 %/K	2,070 %/K
HOSPAL	Hospal	2,000 %/K	2,070 %/K
NIK ALL	Nikkiso Overall-Conductivity	2,050 %/K	2,050 %/K
NIK BIC	Nikkiso Bicarbonate Conductivity	2,200 %/K	2,200 %/K
STANDARD		2,000 %/K	2,070 %/K
USER TC		Adjustable 0 ... 5,000 %/K	Adjustable 0 ... 5,000 %/K

The calculation below shows the drastic effects of an incorrect temperature coefficient.

Example calculation for an incorrect temperature coefficient, using a dialysate with a temperature coefficient of 2.07 %/°C:

Conductivity of solution mS/cm	Temperature of solution °C	Temperature coefficient of meter %/K	Meter display mS/cm	Difference in values %
14,00	37,0	2,07	14,00	0,00
14,00	37,0	1,97	14,17	1,21

### Which temperature coefficient you should use

First of all please double-check the Temperature coefficient with the machine manual.

If you have machines from one manufacturer only in your use the temperature coefficient that the dialysis machine uses for compensation.

If you have different types of dialysis machines in your unit the best solution is to use a temperature coefficient of 2.07 %/°C for **all** machines. This avoids confusion with different readings of the conductivity on different machines.

## Selecting the temperature coefficient



Be aware that changing the temperature coefficient will lead to changes in values displayed by the meter. Therefore only change the temperature coefficient, if you are fully aware of its consequences. If the temperature coefficient setting is wrong it will lead to false display values.

The HDM97 *Pocket* is able to apply dynamic temperature coefficients. It supports a set of 2 coefficients for 25°C and 37°C respective. If only one temperature coefficient is needed it is necessary to set both, the coefficient for 25°C and for 37° to the same value.

## Displaying the temperature coefficient set

To show the current temperature coefficient set press the function key **F** while in conductivity mode. The temperature coefficients are shown with their respective temperatures alternating every 3 seconds. The upper line shows the name of the temperature coefficient set, the middle line shows the temperature and the lower line shows the temperature coefficient associated with the temperature above. If a single temperature coefficient is used the values in the lower line are the same.

The HDM97 *Pocket* enables the user easy to choose between vendor specific temperature coefficients as well as a free adjustable value called "USER TC". To achieve this press the **F** key while in Conductivity / Temperature measuring mode. Use the  and  navigation key to select between the machine manufacturers. By pressing the navigation key **OK** the setting will be applied. Pressing the **ESC** navigation key discards the setting and the meter returns back to measuring mode.

## Adjustment of customer specific temperature coefficient "USER TC"

Pressing the navigation key **OK** while adjusting the temperature coefficient to "USER TC" the HDM97 *Pocket* enters a second dialog to adjust the value itself instead of directly going back to measuring mode.

 → **USER TC**  
**OK** → enters the dialog to adjust the user temperature coefficient set  
**ESC** → cancel and discard setting

Display: **EDIT TC1**  
**37.00 °C** Temperature associated with the coefficient  
**2.071 %/K** (Example 2.070, navigation key  has been pressed once)

Keys:  → increments current value  
 → decrements current value  
**OK** → Applies the value and leads to the adjustment dialog for the second pair  
**ESC** → Discard adjustment and goes on using the last active temperature coefficient setting

Display:	<b>EDIT TC2</b>		
	<b>25.00 °C</b>		Temperature associated with the coefficient
	<b>2.001 %/K</b>		(Example 2.000, navigation key  has been pressed once)
Keys:		→	increments current value
		→	decrements current value
		→	Applies the coefficient set. The displays shows "ADJ OK" for 2 seconds
		→	Discard adjustment and goes on using the last active temperature coefficient setting

### Temperature measurement

This mode is only available if enabled in the basic settings menu. This mode exists to save power in case conductivity measurement is not needed.

By pressing the key  you can switch between unit °C and °F.

### Pressure measurement

In this mode the instrument displays the pressure in the middle line.

#### Unit

By pressing the key  you can choose between units mmHg, mbar, kPa and PSI. The values will be recalculated automatically. It is not necessary to tare or adjust the meter if the unit is changed.

### Offset balancing (tare)

Caused by drift of the pressure probe, the pressure display may sway slightly around zero. The HDM97 *Pocket* is equipped with a tare function which sets the value to zero if drift occurs. While performing this function there must not be any pressure on the pressure transducers inlet. This is assured when the pressure tube is not connected. Then the environmental pressure represents zero.

#### Procedure

1. Make sure there is no pressure applied to the pressure transducer
2. Press and hold down the key  until the upper display line shows "ZERO", this will take three seconds
3. The measuring value is not set to zero.



## Stopwatch (TIMER)

If enabled in the basic settings menu this works as a stopwatch. By pressing the key  e.g. **OK** the stopwatch is being started, by pressing it again it is being stopped. Pressing the key a third time resets the shown time to 00.00.00 and the stopwatch can be used again.



→

**TIMER** die display changes to MM/SS/DS automatically after 2 seconds

**MM/SS/DS**

**00.00.0** ← tenth seconds

↑ ↑

↑ seconds

minutes

After 59 minutes and 59.9 seconds the display form changes so that hours, minutes and seconds are being displayed. The tenth second display is not available then. After 9 hours, 59 minutes and 59 seconds the stopwatch stops itself.

**HH/MM/SS**

**0.00.00** ← seconds

↑ ↑

↑ minutes

hours

## Flow measurement

This mode is available only if a flow sensor is connected.



**Sensor Out**

**Filter Sensor In**

The flow sensor must be connected with the interface shown in the picture below.



Please make sure to connect fluid correctly the sensor. The sensor has a label showing the flow direction.

The head line of the display shows the measuring unit *ml/min* and after a space the unit for the accumulated value as *l* for liter

In the middle line the flow in ml/min and in the bottom line the accumulated value in liter is shown.

For high accuracy measurement we recommend to rinse the flow sensor with 100 ml/min RO water for 15 min.

### Reset accumulated value

The accumulated flow value can be resettled at any time following the procedure below.

Procedure

1. Press and hold down the key  until the upper display line shows "ZERO", this will take three seconds
3. The accumulated value is not set to zero.

### Attenuation of reading

The reading of the current can be attenuated.

With button  the selection menu can be called.

Values up to 180 can be selected. The selected value represents the time in seconds, during the values are averaged.

# Adjustment and Calibration

## Precautions



**Use the adjustment functions only if you are fully aware of their consequences. Wrong adjustment values will lead to false measurement values and in consequence these false values lead to hazards for patients.**

## Definitions

### Adjustment

An adjustment is a process in which properties of a measuring device are changed in that way that its scale or display reads the same as an reference device or standard.

### Calibration

Calibration is the process of establishing the relationship between a measuring device and the units of measure. This is done by comparing a device or the output of an instrument to a standard having known measurement characteristics. For example the length of a stick can be calibrated by comparing it to a standard that has a known length. Once the relationship of the stick to the standard is known the stick can be used to measure the length of other things. Adjustments are checked by performing a calibration.

## General

According to ISO 10012-1 "General Quality Assurance Requirements for Measuring Equipment" the error attributable to calibration should be as small as possible. In most areas of measurement, it should be not more than one third and preferably one tenth of the permissible error of the confirmed equipment when in use. In other words the accuracy of the references to adjust or calibrate the HDM must have at least a tree times higher accuracy.

After a channel has been adjusted it is necessary to calibrate it afterwards. If inconclusive or incomprehensible values are shown in the display it is highly recommended to do a calibration.

The procedure of calibration and adjustment is slightly different for each channel. Proceed with an adjustment only if you have fully understood the procedure.

- Conductivity is adjusted only at one sampling point. No further points are needed due to the high linearity of the conductivity probe.
- The temperature sensor does not drift. Normally there is no adjustment necessary. The adjustment is done by using two sampling points.
- The pressure transducers drift is linear within its tolerances, so there is a tare function (see chapter "Measuring"). Normally there is no adjustment necessary. The adjustment is done by using seven sampling points.
- Due to the kind of their technology it is necessary to adjust pH probes periodically. This is done using two sampling points.

If you are uncertain of entering the correct adjustment values you can discard the values and cancel the adjustment procedure at any time.

## Handling of Reference Solutions

IBP Medical standard solutions are produced under ISO13485:2003 quality management. They are traceable to NIST and PTB Standards Reference Materials and are sealed with tamper-evident packaging.

To ensure standard solution and calibration/verification accuracy

- Keep solutions tightly capped to avoid evaporation
- Do not return used solutions to the storage bottle
- Do not remove solutions from their original bottle
- Keep the solutions in a cool place
- Use only fresh reference solutions for calibration and verification
- Use the solution immediately after pouring, evaporation will cause errors
- Discard solution the appropriate number of days after opening the bottle
- Discard solution after the expiration date

## Adjusting the HDM97 *Pocket*



**Use these functions only if you are fully aware of their consequences. Wrong adjustment values will lead to false measurement values and in consequence these false values may lead to hazards for patients.**

The entire procedure is guided by using software, it is not necessary to open the instrument. To invoke the adjustment menu first select the channel you want to adjust. Then press and hold down the keys  and  at the same time for at least 5 seconds. If you want to cancel the operation press navigation key .

### Temperature

Temperature adjustment is done using 2 sampling points. The order of the adjustment points is preassigned and the value range of each point is restricted by boundaries.

Point	Display	Temperature	Boundaries
1.	"TP1"	25.000 °C	±1.500 °C
2.	"TP2"	55.000 °C	±1.500 °C

Normally an adjustment is not necessary because the sensor has no drift.

The temperature sensor must be dipped into a liquid with known temperature. This known temperature must be adjusted using the navigation keys  and . It is mandatory to wait until the measuring value becomes stable. After the measuring value finally becomes stable the adjustment point can be applied using the navigation key . The instrument jumps immediately to the next adjustment point, so the procedure can be repeated. After the last adjustment point has been applied the instrument displays "CAL OK" in the upper line of its display. If the adjustment is not correct, the instrument displays "CAL ERR" in the display along with a long warning beep. In this case the adjustment values will be discarded and the procedure has to be started from the beginning. A failed adjustment can also be an outcome of a defective instrument.

## Pressure

The pressure transducer has no nonlinear drift so an adjustment is normally not necessary. Use the tare function to balance the pressure transducers environmental pressure change drift.

The pressure channel is adjusted using 7 sampling points. This improves the accuracy of the pressure measurement, because it recalculates slightly nonlinearities of the pressure transducer. The order of the adjustment points starts with point 1 and is preassigned. The value range of each point is restricted by boundaries.

Apply the pressure values to the instruments pressure transducer inlet as shown in the table below. Adjust the value of your reference or standard to the HDM using the navigation keys  and . It is mandatory to wait until the measuring value becomes stable. After the measuring value finally becomes stable the adjustment point can be applied using the navigation key **OK**. The instrument jumps immediately to the next adjustment point, so the procedure can be repeated. After the last adjustment point has been applied the instrument displays "CAL OK" in the upper line of its display. If the adjustment is not correct, the instrument displays "CAL ERR" in the display along with a long warning beep. In this case the adjustment values will be discarded and the procedure has to be started from the beginning. A failed adjustment can also be an outcome of a defective instrument.

Point	Display	Pressure	Boundaries
1.	"PR1"	-500 mmHg	±10 mmHg
2.	"PR2"	-250 mmHg	±10 mmHg
3.	"PR3"	0 mmHg	±10 mmHg
4.	"PR4"	150 mmHg	±10 mmHg
5.	"PR5"	300 mmHg	±10 mmHg
6.	"PR6"	900 mmHg	±10 mmHg
7.	"PR7"	1500 mmHg	±10 mmHg

## Conductivity

The adjustment of the conductivity channel is done by adjusting the cell parameter. To achieve this only one adjustment point is necessary.

Point	Display	Cell parameter	Initial Conductivity
1.	"CD xx.xx°"	0.xxxx	14.000 mS/cm

First of all set the displayed adjustment value to the value of your reference solution. Then make sure that the solution has a stable temperature of exactly 25.00 °C. Wait until both the temperature and the conductivity values in the display become stable before pressing the navigation key **OK**. After that the cell parameter is being calculated and will be stored in the instrument.

## pH

It is necessary to adjust the pH channel before each usage. It is achieved using two adjustment points.

Point	Display	Initial pH value
1.	"PH1"	6.00
2.	"PH2"	8.00

A pH adjustment has to be done according to the following instructions:

Set the mode of temperature compensation to the mode you have to use afterwards. Prepare 2 buffer solutions which should not be more different than 2 pH values from each other and which are around the value you expect from your measurement. If you expect to measure values around pH 6 then use the pH 5 and pH 7 buffer solutions for instance. The adjustment has to begin with the lower pH value and to end with the higher pH value. Preset the buffer solutions value using the ▲ and ▼ navigation keys.

Before dipping the probe into the buffer solution it has to be flushed with distilled water and dried by carefully dabbing it. Depending on the mode you have chosen the temperature sensor must also be flushed and dried. Then dip both into the buffer solution and wait until the measuring values become stable. In this case apply the value using the navigation key **OK**. Flush and dry the probes before continuing with the second adjustment point. The second point is adjusted in the same way. After the last adjustment point has been completed the instrument displays "CAL OK" in the upper line of its display. If the adjustment is not correct, the instrument displays "CAL ERR" in the display along with a long warning beep. In this case the adjustment values will be discarded and the procedure has to be started from the beginning. A failed adjustment can also be an outcome of a defective instrument.

### Cancelling the adjustment procedure

If you are uncertain to have entered the correct values at any step in the adjustment procedure you can cancel the process at any time. Cancelling the operation will discard the adjustment values.

### Restoring adjustment values with factory adjustment values

If you have failed to adjust the instrument properly, you can overwrite the user adjustments with a factory adjustment value set which has been stored into the instrument during production process. Please refer to chapter "Basic settings".

## Maintenance of HDM97 *Pocket*

Generally the HDM97 *Pocket* can be considered as an easy-care instrument. Like all measuring instruments there is a minimum of maintenance and care necessary to ensure that all functions work flawlessly.

Operate the instrument only in a dry environment, and do not touch it with damp hands. Ensure that no fluids intrude into the interior of the device, or into the sockets at the front.

There are no parts in the HDM97 *Pocket* which you can repair yourself. Contact our technical support team in the event of any malfunction. If the device should become damaged, or malfunction, send it to IBP or your distributor for repair.

### **Storage**

Keep the device in a dry place. Suitable is, for example, the original packaging in which you have received the HDM97 *Pocket*, or the carrying case offered by IBP as accessory. If you do not use the device over a longer period of time, you should connect it to the included charging device about every month for one hour to avoid exhausting the batteries.

### **Cleaning**

Never clean the device with any fluids! In case of pollution you can wipe the surface of the HDM97 *Pocket* with a dry and clean cloth. For measurements in the lower conductivity range, the cleaning of the conductivity/temperature measuring cell is recommended. Remove water-soluble substances by rinsing with deionized water, fats and oils with warm water and household dish washing detergent. Lime and hydroxid crusts can be dissolved by a 10% citric acid solution. In each case, the measuring cell needs to be washed with deionized water after the cleaning. Basically, the conductivity measuring cell does not decay over time. Particular measuring media (for example, strong acids and caustic solutions, organic solvents), or too high temperatures shorten the service life considerably, and/or lead to damage.

### **Caring for the conductivity/temperature electrode**

Thorough cleaning of the electrode is particularly important for measuring low conductivities. Water-soluble substances must be removed by rinsing with deionized water. Remove lime deposits with a 10% citric acid solution. To do this, immerse the electrode in the solution for 2 hours at room temperature, then rinse it thoroughly with deionized water. Do not touch the electrode surface with your fingers. If necessary, clean the surface with acetone.

### **Caring for the flow sensor**

After use make sure to rinse the flow sensor with deionized water for a couple of minutes. Please clean the inlet filter on a regular basis.

## Calibration period

When delivered new from the factory, each device is adjusted and calibrated before shipment. To ensure reliability, we recommend the following intervals for verifying calibration:

<b>Parameter</b>	<b>Calibration period</b>
Temperature	Once per year
Conductivity	Once per year
Pressure	Once per year
Flow	Once per year
pH	Before each use

***A complete function test and calibration of the device should be done once a year.***

According to ISO 10012-1 "General Quality Assurance Requirements for Measuring Equipment" the error attributable to calibration should be as small as possible. In most areas of measurement, it should be not more than one third and preferably one tenth of the permissible error of the confirmed equipment when in use. In other words the accuracy of the references to adjust or calibrate the HDM must have at least a three times higher accuracy.

***For liability reasons you should not carry out these tests and calibration and verification yourself. If IBP performs the calibration, you will receive a calibration certificate in accordance to ISO9001 which documents the calibration/verification results.***

## Tolerances

For calibration, the instrument must meet the deviations listed below.

Channel	Reference Value	Allowed deviation		Notes
Temperature	25,00 °C	± 0,05 °C		
	34,00 °C	± 0,05 °C		
	37,00 °C	± 0,05 °C		
	40,00 °C	± 0,05 °C		
	55,00 °C	± 0,07 °C		
	80,00 °C	± 0,07 °C		
Conductivity	74,00 µS/cm	± 0,60 µS/cm		
	147 µS/cm	± 0,60 µS/cm		
	720 µS/cm	± 6,00 µS/cm		
	1410 µS/cm	± 6,00 µS/cm		
	2,77 mS/cm	± 0,03 mS/cm		
	6,70 mS/cm	± 0,03 mS/cm		
	12,88 mS/cm	± 0,03 mS/cm		
	16,00 mS/cm	± 0,03 mS/cm		
Pressure		General	BM & Opt.H	
	+ 1900,00 mmHg	± 2 mmHg	± 2 mmHg	
	+ 1500,00 mmHg	± 2 mmHg	± 2 mmHg	
	+ 1000,0 mmHg	± 2 mmHg	± 1 mmHg	
	+ 300,0 mmHg	± 1 mmHg	± 0,5 mmHg	
	+ 200,0 mmHg	± 1 mmHg	± 0,5 mmHg	
	+ 100,0 mmHg	± 1 mmHg	± 0,5 mmHg	
	0,0 mmHg	± 1 mmHg	± 0,5 mmHg	Open inlet
	- 100,0 mmHg	± 2 mmHg	± 2 mmHg	
	- 400,0 mmHg	± 2 mmHg	± 2 mmHg	
- 600,00 mmHg	± 2 mmHg	± 2 mmHg		
pH	355,2 mV = pH 1	± 0,02		$\Delta E = -59,2 \text{ mV} / \Delta \text{pH}$
	177,6 mV = pH 4	± 0,02		$\Delta E = -59,2 \text{ mV} / \Delta \text{pH}$
	118,4 mV = pH 5	± 0,02		$\Delta E = -59,2 \text{ mV} / \Delta \text{pH}$
	59,2 mV = pH 6	± 0,02		$\Delta E = -59,2 \text{ mV} / \Delta \text{pH}$
	0 mV = pH 7	± 0,02		$\Delta E = -59,2 \text{ mV} / \Delta \text{pH}$
	- 59,2 mV = pH 8	± 0,02		$\Delta E = -59,2 \text{ mV} / \Delta \text{pH}$
	-118,4 mV = pH 9	± 0,02		$\Delta E = -59,2 \text{ mV} / \Delta \text{pH}$
	-177,6 mV = pH 10	± 0,02		$\Delta E = -59,2 \text{ mV} / \Delta \text{pH}$
	-355,2 mV = pH 14	± 0,02		$\Delta E = -59,2 \text{ mV} / \Delta \text{pH}$

## Troubleshooting

**Problem:**

USB cable connected, but the battery is not being charged

**Reason:** USB port of laptop does not provide enough power

**Solution:** Connect the meter to the battery charger or a USB port which is able to supply at least 500 mA.

**Reason:** USB cable broken

**Solution:** Replace cable

---

**Problem:**

The meter can not be switched on

**Reason:** Battery is empty.

**Solution:** Connect the meter to a battery charger or an USB port.

**Reason:** Other problems.

**Solution:** Press and hold the on/off key for at least 5 seconds. The meter will perform a "Low-Level-Reset" and should start afterwards.

---

**Problem:**

The meter powers off itself after a couple of seconds.

**Reason:** Battery empty.

**Solution:** Connect the meter to a battery charger or an USB port.

---

**Problem:**

pH channel can not be selected (only Version *BP*)

**Reason:** pH channel is disabled in basic settings menu.

**Solution:** enable pH channel in basic settings menu.

---

**Problem:**

Pressure value does not tare / zero balancing does not work (only versions *BO* and *BP*)

**Reason:** There is still pressure applied to the pressure transducer.

**Solution:** Disconnect pressure hose directly at meter.

## Technical data

### Conductivity

0 to 30,0 mS/cm, 4-Pole probe,  
Resolution and accuracy:

Range	Resolution	Accuracy
0.0 – 200 $\mu$ S/cm	0,1 $\mu$ S/cm	$\pm 0,6 \mu$ S/cm
201 - 1999 $\mu$ S/cm	1 $\mu$ S/cm	$\pm 6\mu$ S/cm
2,00 – 30 mS/cm	0,01 mS/cm	$\pm 0,03$ mS/cm

Reference temperature 25° C

Temperature compensation is 2,070 % per Kelvin  
Adjustable from 0.00 to 4,00 %/Kelvin  
Or selectable from dialysis machine manufacturer preset values.

### Temperature (only versions BN, BO, BP and BQ)

Range 0 ... 100 °C,  
Resolution 0,01 °C  
Accuracy 25 ... 40°C  $\pm 0,05^\circ\text{C}$  , otherwise  $\pm 0,1^\circ\text{C}$

### Pressure (BO, BP and BQ versions, fluid resistant ceramic sensor)

Units mmHg, kPa, mbar, PSI  
Range -700 ... 1900 mmHg, 2.5 Bar  
Resolution 0,1 mmHg  
Accuracy 0 ... 300 mmHg  $\pm 1$  mmHg  
Otherwise  $\pm 2$  mmHg  
Fluid resistance yes

### Pressure (only Version BM and Option H for BQ Version)

Units mmHg, kPa, mbar, PSI  
Range -700 ... 1900 mmHg, 2.5 Bar  
Resolution 0.01 mmHg  
Accuracy 0 ... 300 mmHg  $\pm 0.5$  mmHg  
301 ... 1000 mmHg  $\pm 1$  mmHg  
Otherwise  $\pm 1$  mmHg  
Fluid resistance no

### pH (only BP Version)

Range pH 0 ... 14 ; max. slew rate 70mV (118%) /  $\Delta$ pH  
Resolution pH 0,01  
Accuracy pH  $\pm 0,02$   
DC input impedance min. 10 Tera Ohm (min.  $10 * 10^{12}$  Ohm)  
Input leakage current typ.  $\pm 2$  fA ; max.  $\pm 2$  pA  
Temperature compensation with temperature measurement or preset values to choose from.

### Flow (only version BL and BQ)

Range 100 ... 2500 ml/min  
Resolution 1 ml/min  
Temperature range: 0 ... 85°C  
Accuracy 100 ... 250 ml/min  $\pm 3\%$  from current value,  
otherwise  $\pm 3\%$  from upper range value

**Internal power source**

Internal Li-Ion-Polymer battery, rechargeable with USB 5V DC / max. 500mA  
Operating time, depending in channel, up to 60 hours.

Nominal Voltage	3,7 V
Max. charge voltage	4,2 V (internally generated an monitored)
Capacity	850 mAh or 3,1 Wh
Battery charge time	approx. 3 hours
Device use time	estimated 40 hrs (displaying conductivity) up to 120 hrs (stopwatch) USB not connected

**External power supply**

Input voltage	100-240 V AC/ 50-60Hz
Output voltage/current	5 V DC, 500 mA min.
Plug	Mini USB A-Plug

**Degree of protection**

Meter	IPX1
Conductivity probe	IP66 , max. 0,5 m immersion depth, max. 5 min

**Size and weight**

Size	165 x 82 x 33 mm
Weight	400 gr. without battery charger, including conductivity/temperature probe

These technical data may be changed without notice!

## PC Interface

### USB Interface

The USB interface of the HDM97 *Pocket* is electrically isolated from the appliance part. Two drivers are needed for installation. One driver is made for invoking the interface like any other USB device and the other driver is used to emulate a serial COM port device.

### Interface parameters for serial COM port emulation

The HDM97 *Pocket* uses the preset values below for serial COM port emulation:


<b>Property</b>	<b>Value</b>
Baud rate	9600 bps
Data bits	8
Parity	None
Stop Bit	1
Handshake	none

Driver installation, example for *Microsoft® Windows® XP Service Pack 2*

1. Insert the Installation-CD in the CD or DVD drive of your computer.
2. Connect the HDM97 *Pocket* via the USB-cable with a free USB port of the running computer. It does not matter whether the instrument is switched on or off.
3. A "Found New Hardware Wizard" Window will appear on the screen like that right here from the text.
4. Choose "No, not this time" and afterward click on the button "Next".




5. Driver source selection dialog: Choose "Install from a list or specific location (Advanced)", and click on the "Next" button afterwards

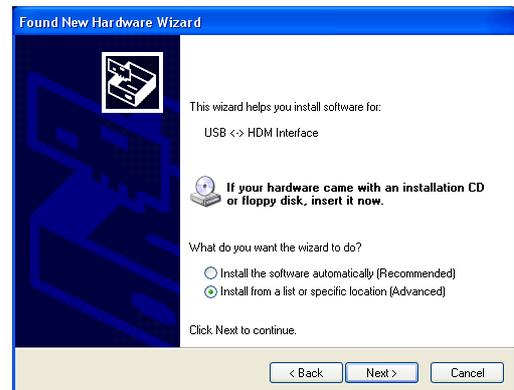

6. Choose "Search for the best driver in these locations" and check "Include this location in the search:". Then use the "Browse" button and choose your CD/DVD drive's letter and add the folder "drivers" to it. In the given example on the right side the CD/DVD driver has drive letter D:\ assigned.


Click on the "Next" button afterwards.

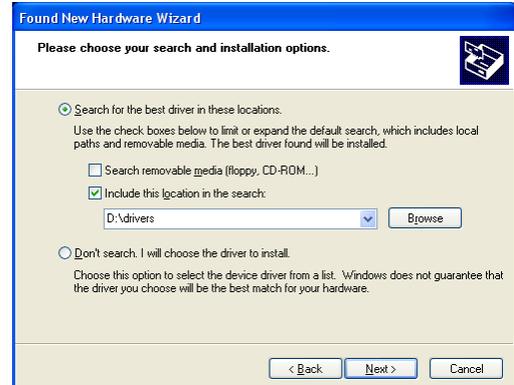

7. Windows Logo Test : Choose "Continue Anyway".



8. The driver installation is in progress. This can take a couple of minutes.

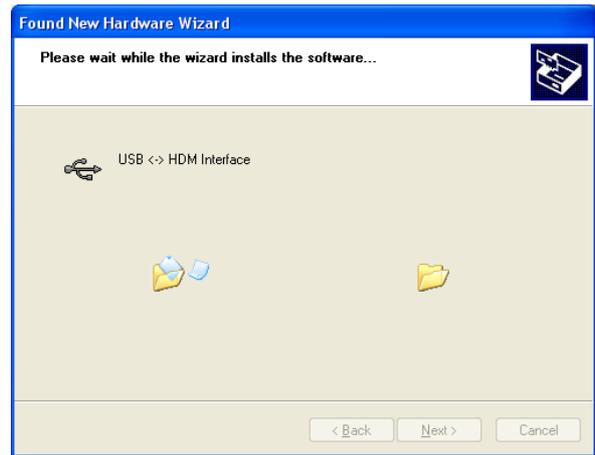

9. The driver installation was successful, click on the button "Finish" to complete the procedure.




After you have finished this procedure another similar Wizard will popup. Do exactly the same you did in the first procedure as described earlier. One installation is for the COM port emulation, the other installation is for the USB interface itself.


## Checking the driver installation (optional)

The USB interface of the HDM97*Pocket* emulates a serial COM port and offers a proprietary interface to communicate directly with the USB interface.

Use the "My Computer/Manage" function to open the Device Manager and search for a Group called "Ports (COM and LPT)". You should find an entry called "USB <-> HDM Interface". At the end of the line you will find the number of the associated COM port (in this case "COM3").

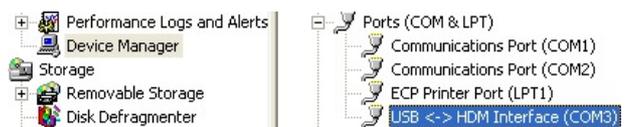

In addition you will find the USB communication interface in the Group "USB-Controller" It is named " USB <-> HDM Interface ".

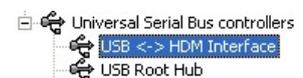

## Data transfer

To receive measurements from the HDM97 *Pocket* via the USB interface, the following abbreviations must be sent via the interface. The entire communication takes place in readable ASCII-characters.

<b>Command to HDM97 <i>Pocket</i></b>	<b>Reaction of the HDM97 <i>Pocket</i></b>
HDMCMO	activates the data transfer
HDMCMF	deactivates the data transfer
HDMEIH	transmits all units
val	transmits the values

## Format of the values

Conductivity in S/cm	9 signs
Temperature in °C	7 signs
Pressure in mmHg	9 signs
pH	5 signs only version <i>BP</i>
Flow in ml/min	5 signs only version <i>BQ</i>
CR, LF	2 signs

The separation of the values is done by using "/"

## Example

Command „hdm“ – HDM97*BP* sends: 0.1234567/123.123/ -12345.12/12.12 + CR LF

Command „HDMEIH“ – HDM97*BP* sends: S/cm,°C,mmHg,pH CR LF

# IBPView

## What is IBPView

**IBPView** can be obtained optionally for the **HDM97Pocket** and can be used as well for other HDM instruments. A demo version is included in the folder \Software\ on the driver CD. This demo can not save, export or print and the acquisition time is limited to 30 seconds.

**IBPView** is a data acquisition software, which records and plots measured values from a **HDM97Pocket** device. It works with a USB connection of the **HDM97Pocket**. The software can also be used with several hardware devices, such multimeters from *Voltcraft* or *Metex*.

**IBPView** is compatible with *Microsoft® Windows® 2000*, *Microsoft® Windows® XP* and *Microsoft® Windows® Vista®*

Several curves can be plotted into a single diagram. Several diagrams can be arranged on one or more pages if necessary.

Plot options like pen width or pen color can be used to create perfectly styled diagrams. You can even create your own plotter designs.

Comfortable zoom functions are available on a single click to analyse plot details or to return to the plots overview.

Setting markers makes it easy to read differential values of a plot, or to determine the amplitude or period of a sampled signal.

A comfortable preview allows you to add comments or project data when printing diagrams.

Export functions can be used to use measured data with other applications like MS-Excel, etc.

The help function of **IBPView** includes instructions for basic operations using a small project as an example. If you follow these instructions it will be a good start for you in understanding the basics.

## Installation

Insert the CD „IBPView“ into the CD-Drive. The installation will start automatically. In case open the directory „SETUP“ and start setup.exe. On then first menu select the language, the installation process is self explaining.


## Quickstart

On the CD „**HDM97Pocket**“ you will find templates for all versions of the **HDM97Pocket** . These templates are prepared for two, three and four measuring channels. Do not use a template which is prepared for more measuring channels than measuring channels meter has.

## Disconnect USB

Before terminating the USB connection please terminate the IBPView software. Disconnecting with running data acquisition may end in a system freeze.

## Main Window

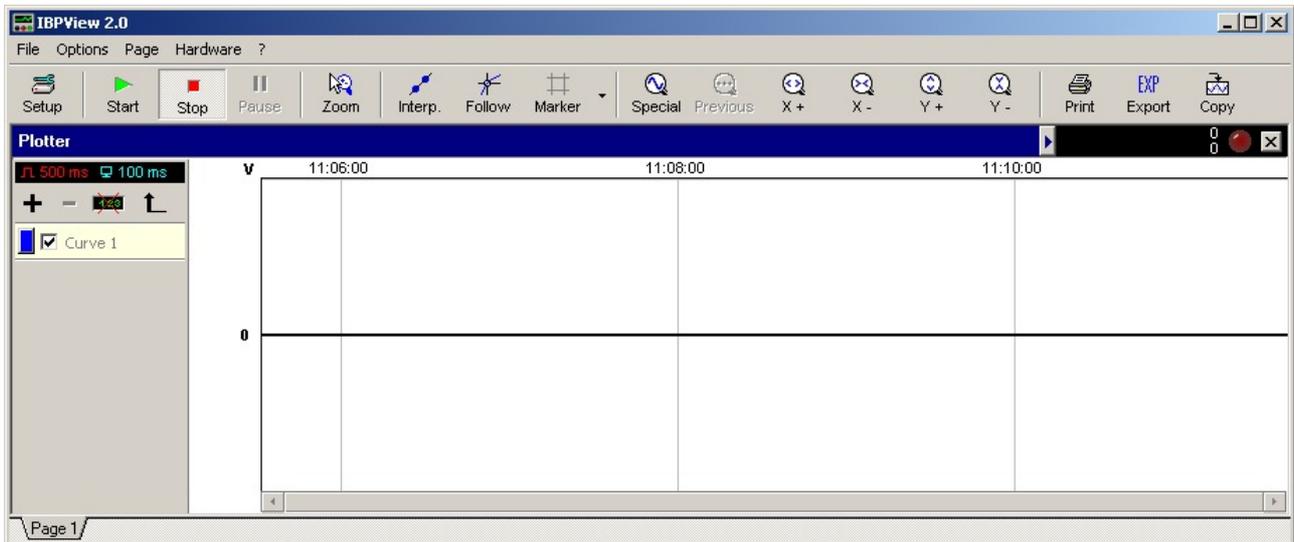

## General

This manual describes only the basic functions which are necessary to get familiar with **IBPView**. Further information can be obtained using the Software's help function. **IBPView** is intentionally not limited to HDM instruments. It can be configured independently and therefore achieves its maximum flexibility to be ready for future IBP products.

## Hardware Interface

**IBPView** uses serial COM ports for communication. This can be common RS232 serial interfaces like many multimeters on the market use, or, like the **HDM97Pocket** virtual serial ports based on USB. Using virtual COM ports it is necessary to define the number of the port once (COM1, COM2, COM3, etc.). The COM port number of the hardware is defined automatically during the driver installation. The application must know this number, in order to communicate with the device. For obtaining the number for the **HDM97Pocket** please refer to chapter "Checking the driver installation". In this example "COM3" is used. Be aware that connecting the **HDM97Pocket** to different physical USB receptacles will result in different COM port numbers, one for each receptacle. This is done to make multiple HDM meters work with one computer concurrently.


## Template projects

Project files are named in the way "[Name].ibp". To get an easy start with **IBPView** there are preconfigured version specific templates for each type of instrument. These templates include all common setting for start working with the application. Only the number of the COM port must be adjusted to the number the driver installation has chosen if it is not coincidentally the same number is used. For help to select the right COM port read also chapter "Checking the driver installation".


Templates are available on the CD "HDM97 Pocket".

## Selecting hardware

1. Click on "Hardware" in the main window and choose menu entry "Define...". Then a window named "Hardware configuration" like that right from this text will open. Open in the left box "Supported hardware devices" the group "IBP", click on the entry "IBP Universal" and then drag and drop it to the right box "Existing hardware devices".

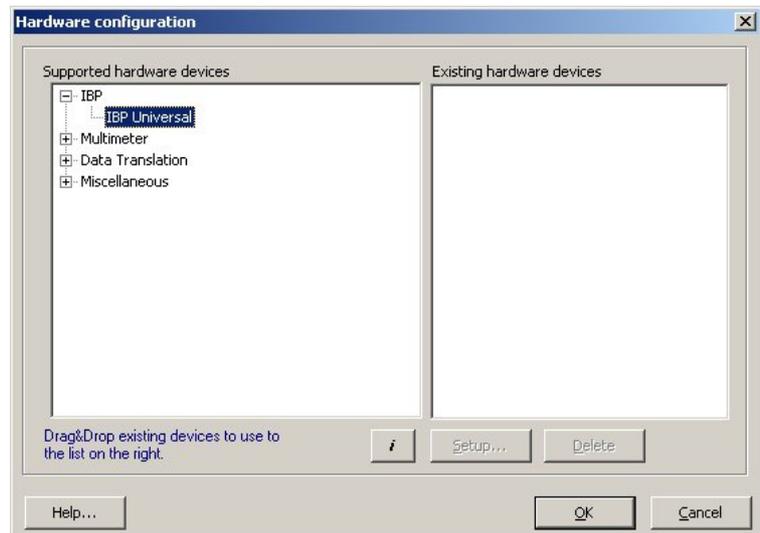


2. After the drag and drop procedure a window opens which can be used to set the number of the serial COM port and the number of channels for the instrument. For a HDM97BP four channels are possible. In case the pH channel is disabled in the base settings of the HDM97BP, the meter has only three measuring channels. Selecting of a wrong channel number than the meter really provides will result in a not working data transmission.

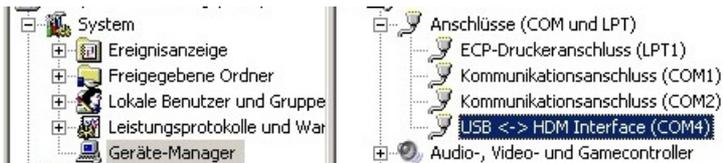


To find out which COM-Port the USB-Driver has selected please select the Windows System Manager and than the Device Manager. Under group "Connctions (COM and LPT)" you should find the port.

3. After this the device is displayed in the right box. Other hardware devices can be added if needed. Click on "OK" to close the window.

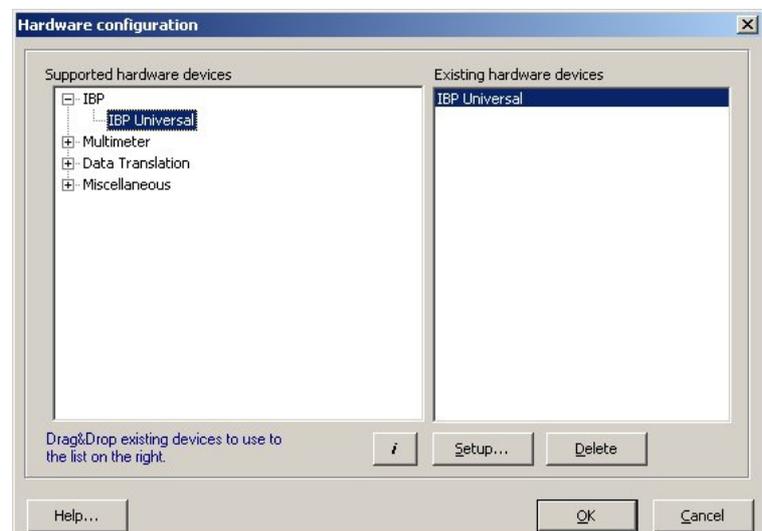


### Assigning graphs to hardware channels

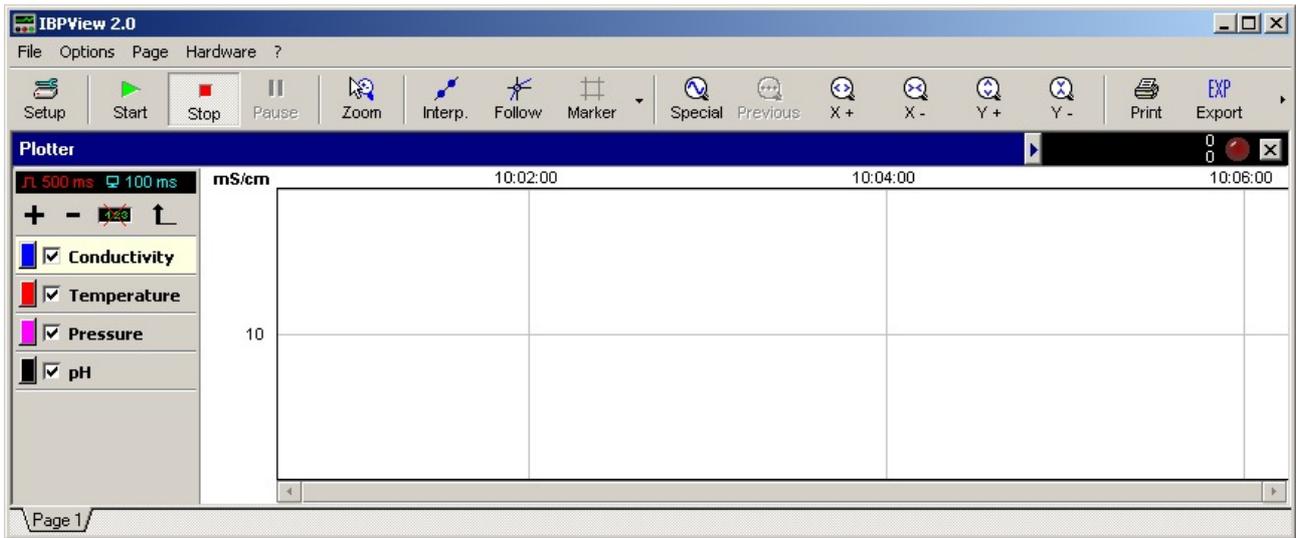

In the main window the list of channels is located left. The assignment of a particular graph to a specific hardware channel is done by performing a double click on the channel's name in the list.

This double click leads to a dialog "Channel configuration" (picture in the middle). You can change Name, Basic units, Decades, Ranges and much more. By clicking on the button "Assign" you can assign each graph a hardware channel, like shown in the picture right.

For conductivity please use the base unit S/cm and select for example "m (milli)" as Decade to show the values in mS/cm.

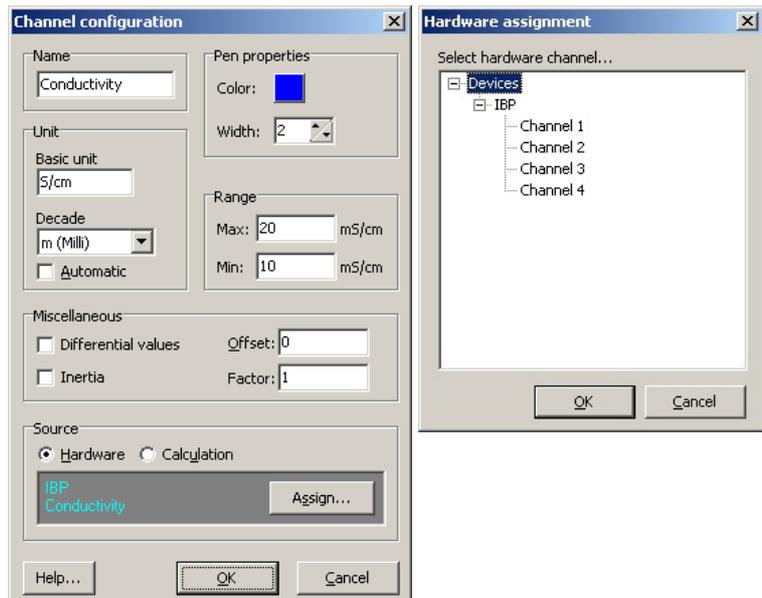

## List of channels

Depending on the version the HDM97 *Pocket* is equipped with up to 4 channels

Ch.	Description	Range	Unit	HDM97 <i>BL</i>	HDM97 <i>BM</i>	HDM97 <i>BN</i>	HDM97 <i>BO</i>	HDM97 <i>BP</i>	HDM97 <i>BQ</i>
1.	Conductivity	0.000 .. 30.00	mS/cm			✓	✓	✓	✓
2.	Temperature	0,00 .. 100,00	°C			✓	✓	✓	✓
3.	Pressure	-750,0 .. 1900	mmHg		✓		✓	✓	✓
4.	pH	0 .. 14.00	without					✓ If not disabled	
4.	Flow	0 ... 5000	ml/min	✓					✓ If not disabled

If the wrong number of channels is selected the data transmission will not work correctly. In case the pH channel is disabled in the base settings of the *HDM97BP*, the meter has only three measuring channels.