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



Switch on the instrument

The power switch is on the right side of the instrument.

Start screen shows: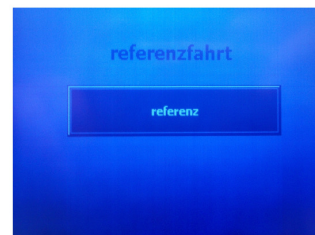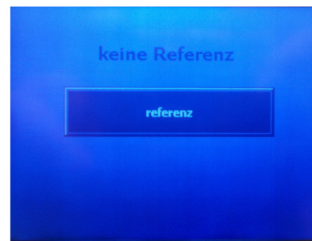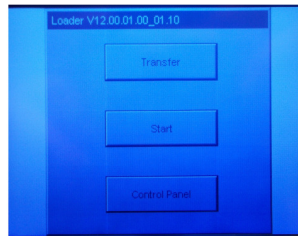

Press reference

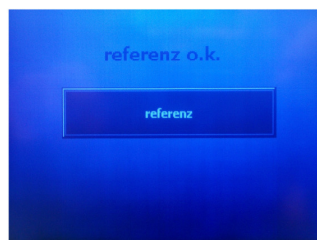
selfcontrol starts

Press reference

Press F2 / Program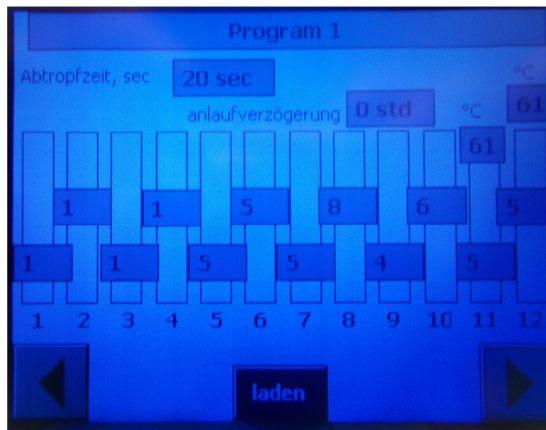

The liquid container have numbers from 1 – 12


Choose the processing time for container 1 by pressing the grey field, insert the processing time for container 1 for example 60 minutes.

← 60 Minutes


Confirm by pressing ENTER ↑

Insert time for the next container, If there is no time / 0 minutes the basket will skip this container. Insert time for all the containers.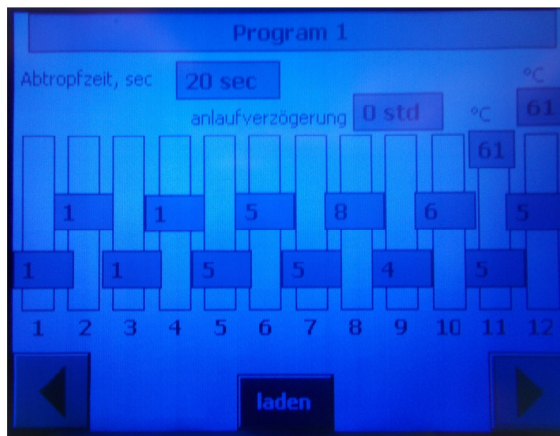

← Choose the temperature
Confirm by pressing
ENTER

If desired insert, the start time delay for weekends or holidays ,
this time is set in hours not in minutes.
The maximum delay is 4 days .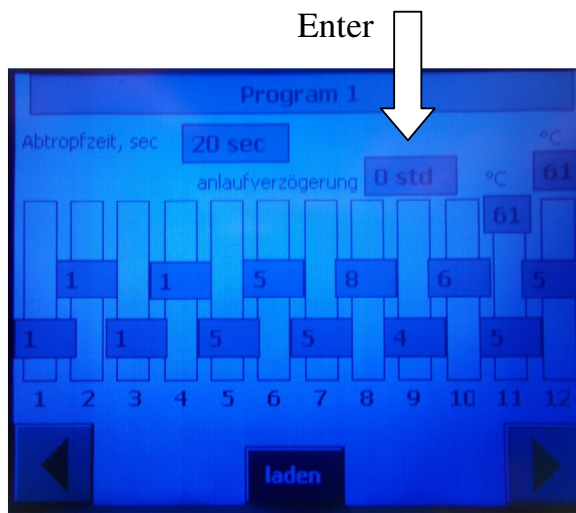

After entering all parameters , press laden (load), the program will
stored .


Fill all containers with the appropriate solvent / paraffin .

The container should be filled up to a maximum 4 cm below the rim .

If the paraffin is not liquid , it may take up to 5 hours until it
is completely melted .

Hang the basket with the tissue cassette into the holder ,
Close the door .

Press F1 / Auto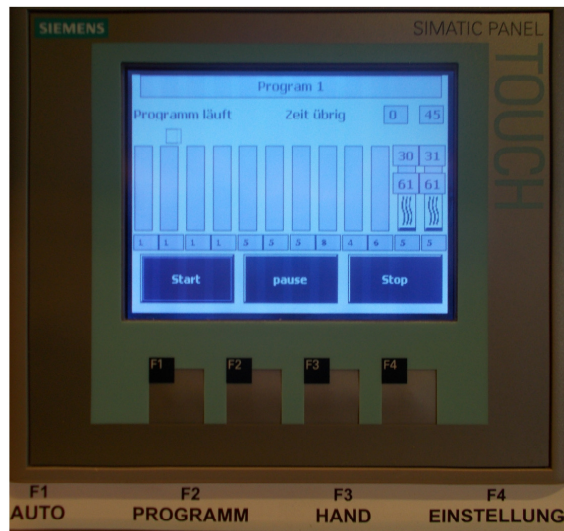


Press Start the program starts..

After pressing start do not open the door again !

If you open the door the basket will return to the start position.

After completion of the program, press F3 / HAND , you can now lift the basket with the arrow keys up .

Take out the basket. Press the Home button , the device moves into the Start position .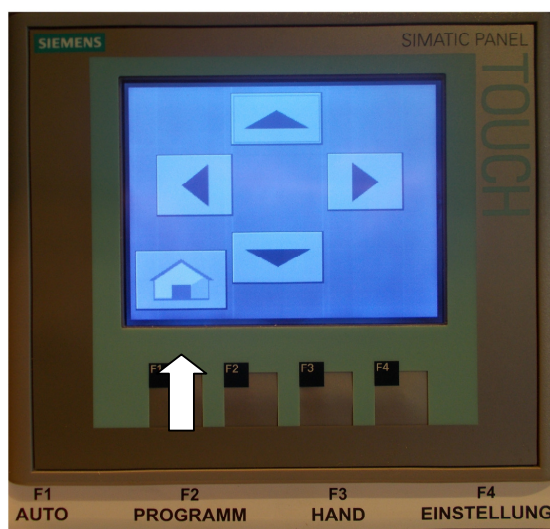

If you want to start a further program , move the lid down by using the arrow down key .

You can save up to 5 programs

If the wax does not reach the intended temperature the basket remains in the last container standing in front of the paraffin .

In case of power failure , the basket always remains in liquid and must be continued manually.

The device has a battery backup which the device enables to continue to operate for up to 2 minutes.