

iFLEX-iRIS™

Laser Systems

The Qioptiq iFLEX-iRIS™ solid state laser systems offer high performance stability with low amplitude noise in a miniaturized package. They are ideally suited for integration into instruments that demand high performance yet need to retain a small form factor. The lasers are mode-hop free and wavelength stabilized as a result of active temperature control.

All CW iFLEX-iRIS lasers operate in automatic power control mode using an internal feedback loop. This feature provides highly stable output power, and ensures the high power stability performance is maintained throughout the product lifetime.

Well controlled digital modulation can be offered through the electrical interface PINs. This feature must be selected with order.

To ensure optimum performance is maintained throughout the product lifetime, the iFLEX-iRIS laser design is robustly tested to Telcordia specifications.


Product features include:

- · High stability operation
- Low noise
- Miniaturized for easy integration
- Integrated electronics and TEC temperature stabilized control
- Conductively cooled through baseplate
- Automatic Power Control
- Mechanical shutter

Options include:

- Digital modulation
- CDRH compliant interlock unit
- Adaptor cable when an iFLEX-iRIS is used as a retrofit replacement for an iFLEX2000™ laser
- Industry compatible footprints and beam exit positions
- OEM versions on request
- Fiber coupling please refer to the fiber coupled datasheet for information about fiber coupled iFLEX-iRIS lasers



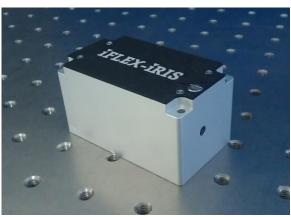



iFLEX-iRIS laser specification overview

Operating Wavelengths

Wavelength (nm)	375	405	445	458	473	488	505	515	520	633	637	642	660	670	730	780	830	852
Laser Power (mW)	20 40 50	50 100 200 220	20 50 75	20 70	75	20 40 100 140	50	20 50 60	30	30 70	20 100	20 40 100	80	10	20	70	100	35

Technical Specifications

Lacor parameters						
Laser parameters						
Center wavelength	± 5*	nm				
Power stability (8 hours)	< 1	%				
Optical Noise (20Hz-2MHz)	< 0.1**	% rms				
Spatial mode	TEM ₀₀	-				
Beam diameter at 1/e ²	0.7 ± 0.2	mm				
Beam divergence (full angle)	Typ < 1**	mrad				
M ² (beam quality)	Typ 1.2	-				
Pointing stability	< 5 µ	ırad/°C				
Polarization ratio	≥ 200:1	-				
Static alignment						
Beam position	< 0.3	mm				
Beam angle	< 5	mrad				
Electrical						
Power supply	12V DC, 1A	-				
Max. base plate temp.	40	°C				
Max. heat dissipation	12	W				
Operation Modes						
CW	0.1 - 100% and true off					
Digital Modulation	TTL signal	-				
Rise time (10% to 90%)	< 1	μsec				
Fall time (90% to 10%)	< 1	μsec				
Modulation bandwidth	DC to 500	kHz				
Laser Dimensions						
Laser only	70(L) x 40(W) x 38(H)	mm				

Environmental conditions					
Storage temperature	10 to 50	°C			
Operating pressure	Atmospheric	-			
Operating temperature	10 to 40	°C			
Operating & storage humidity	Non-condensing, dew point <23°C				

^{*}Center wavelength exceptions at: 637nm (+6/-5nm), 642nm (+5/-7nm), 660nm (+4/-6nm), 670nm (\pm 10nm), and 730nm (\pm 10nm)

^{**}Model specific please contact Qioptiq for details Custom OEM versions on request



iFLEX-iRIS laser specification overview

Operating Wavelengths:

Wavelength (nm)	532	561	594
Laser Power (mW)	20, 40	20, 50	20

Technical Specifications

_							
Laser parameters							
Center wavelength	± 2	nm					
Power stability (8 hours)	< 2	%					
Optical Noise (20Hz-2MHz)	< 0.3*	% rms					
Spatial mode	TEM ₀₀	-					
Beam diameter at 1/e ²	0.7 ± 0.2	mm					
Beam divergence (full angle)	Typ < 1*	mrad					
M ² (beam quality)	Typ ≤ 1.2	-					
Pointing stability	< 5	µrad/°C					
Polarization ratio	> 200:1	-					
Static alignment							
Beam position	< ± 0.3	mm					
Beam angle	< 5	mrad					
Electrical							
Power supply	12V DC, 1A	-					
Max. base plate temp.	40	°C					
Max. heat dissipation	12	W					
Operation Modes	Operation Modes						
CW	50 - 100% adjustment and true off						
Laser Dimensions							
Laser only	70(L) x 40(W) x 38(H) mm					

Environmental conditions					
Storage temp.	10 to 50	°C			
Operating pressure	Atmospheric	-			
Operating temp.	10 to 40	°C			
Operating & storage humidity	Non-condensing, dew point <23°C				


^{*}Model specific please contact Qioptiq for details Custom OEM versions on request



Model: iFLEX-iRIS-X0

iFLEX-iRIS laser with central laser aperture position


Note: it is possible to attach a kineFLEX fiber delivery system to the iFLEX-iRIS-X0 laser with central laser aperture (beam exit) position. The kineFLEX can be attached by customers, as a future upgrade option.

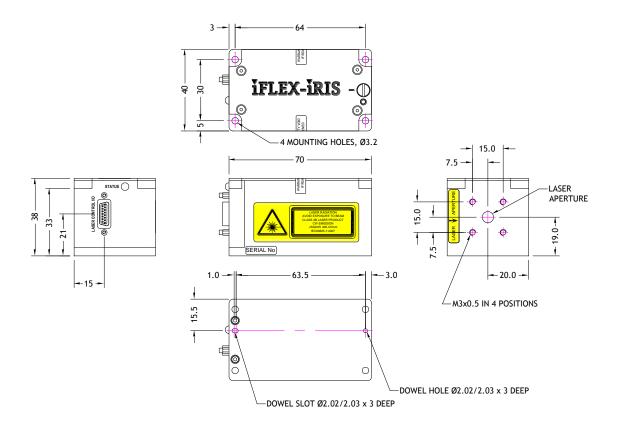

iFLEX-iRIS electrical interface

Pin Function

- 1 Monitor Photodiode Output (Vmon)
- 2 Operating Current Output (Vop)
- 3 External Power Control
- 4 Laser Enable
- 5 Temperature OK Signal
- 6 Reserved
- 7 Laser Supply Voltage
- 8 Reserved
- 9 Reserved
- 10 Reserved
- 11 Chassis Ground
- 12 Digital Modulation Input
- 13 Reserved
- 14 OV Laser Supply
- 15 OV Signal Return

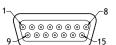


Please refer to iFLEX-iRIS user manual for detailed instructions on electrical control via PINs.



Model: iFLEX-iRIS-X2

iFLEX-iRIS laser with offset laser aperture position

Note: the kineFLEX fiber delivery system cannot be directly attached to the iFLEX-iRIS-X2 laser with offset laser aperture (beam exit) position, and an adaptor interface would be required.

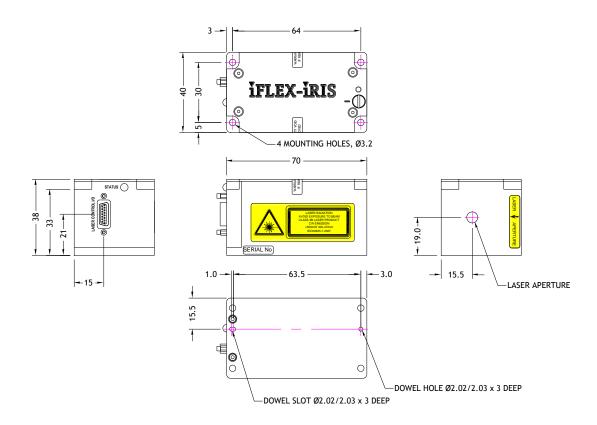


Other industry footprints and beam offsets may be available. Please contact Qioptiq Sales for details.

iFLEX-iRIS electrical interface

Pin Function

- 1 Monitor Photodiode Output (Vmon)
- 2 Operating Current Output (Vop)
- 3 External Power Control
- 4 Laser Enable
- 5 Temperature OK Signal
- 6 Reserved
- 7 Laser Supply Voltage
- 8 Reserved
- 9 Reserved
- 10 Reserved
- 11 Chassis Ground
- 12 Digital Modulation Input
- 13 Reserved
- 14 OV Laser Supply
- 15 OV Signal Return

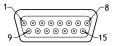


Please refer to iFLEX-iRIS user manual for detailed instructions on electrical control via PINs.




iFLEX-iRIS CDRH compliant interlock remote unit


For CDRH compliance a bench top standalone laser system must be used with a CDRH compliant interlock remote unit power supply. Qioptiq is pleased to offer this iFLEX-iRIS CDRH compliant interlock remote unit for use with the iFLEX-iRIS laser systems.

Note: The iFLEX-iRIS CDRH compliant interlock remote unit includes keyswitch, keys, interlock plug and power supply. This CDRH compliant interlock unit is recommended for bench top use.

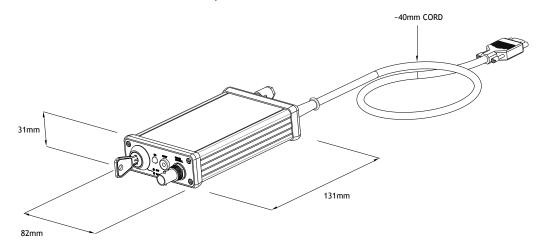


End panel views of the iFLEX-iRIS™ interlock remote unit detailing the location of the different connectors and indicators

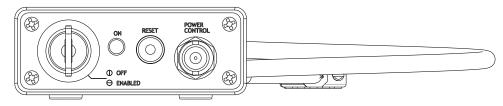

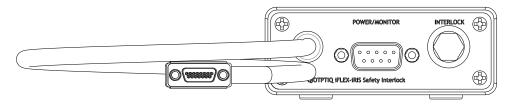

iFLEX-iRIS OEM power supply

When iFLEX-iRIS lasers are integrated into OEM instrumentation then power may be supplied directly from the instrument or from an OEM power supply. Qioptiq is pleased to offer this OEM power supply for the iFLEX-iRIS laser systems.

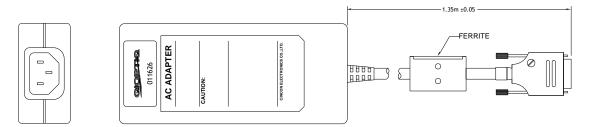


iFLEX-iRIS Adaptor Cable

This allows an iFLEX-iRIS to be installed in place of an iFLEX2000, as a retrofit laser system.



iFLEX-iRIS Laser Order Codes

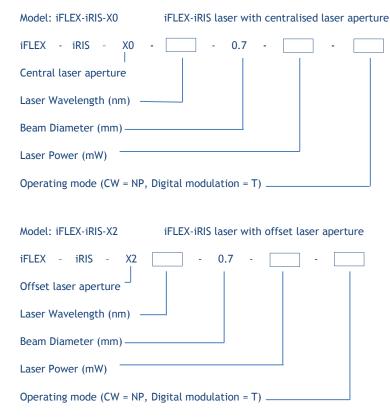

For example: iFLEX-iRIS-X2-640-0.7-50-T

This iFLEX-iRIS laser operates at 640nm and outputs 50mW power. The output beam is collimated to 0.7mm diameter. The beam exit position is offset. This laser may be modulated in current control mode up to 500kHz. This does not include a power supply.

Please remember to select the appropriate power supply for your application

iFLEX-iRIS Power Supply Order Codes

Model: iFLEX-iRIS-CDRH-PSU-SET iFLEX-iRIS CDRH interlock remote unit with power supply, keys and interlock plug

Model: iFLEX-iRIS-OEM-PSU iFLEX-iRIS OEM power supply

Model: iFLEX-iRIS-ADAPTOR-2000 Adaptor cable from iFLEX-iRIS laser to iFLEX-2000 power supply

Note: IEC power lead to wall socket is country specific and may be sourced locally through most electronic retailers. It is not included.

These specifications may be amended without notice under our program of continual improvement, and Qioptiq reserve the right to amend these specifications without notice.

These lasers are designed and manufactured in a factory with ISO9001:2008 certification. These lasers are CE certified. These lasers meet the requirements for IEC60825-1. Only lasers used with a CDRH compliant interlock unit meet CDRH requirements. The addition of a fiber optic output can increase laser hazard. Please follow recommended laser safety guidelines when using these lasers.

For technical information contact:

Qioptiq Photonics Ltd. sales@qpl.qioptiq.com phone +44 (0) 2380 744 500 www.qioptiq.com







iFLEX-iRIS™ and iFLEX-2000™ are trademarks of Qioptiq Photonics Ltd. Copyright ©2014 Qioptiq Photonics Ltd.

