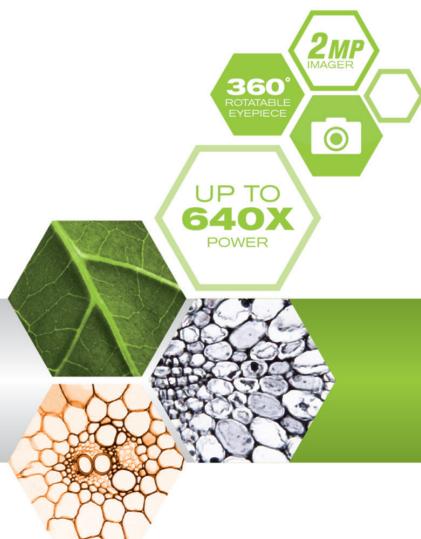
MICRO 360™ +

Multi-Purpose Microscope with 2MP Imager

INSTRUCTION MANUAL MODEL #44126

USB

INTRODUCTION

Congratulations on your Celestron microscope purchase. Your new Micro360+ Dual Purpose Microscope with 2MP Imager is a precision optical instrument made of high quality materials to ensure durability and long life. It is designed to give you a lifetime of enjoyment with minimal maintenance.

Before attempting to use your Micro360+, please read through these instructions thoroughly to familiarize yourself with its functions and operations to maximize your enjoyment and usage. Refer to the microscope diagram to locate parts discussed in this manual.

Micro360+ provides powers from 40x to 640x. It is ideally suited for examining specimen slides of yeasts and molds, cultures, plant and animal parts, fibers,

bacteria, etc. The included 2MP imager is made of high quality materials and can be used with virtually any traditional microscope with 23 mm or 30 mm eyepiece tubes. Please read through the complete instructions to familiarize yourself with all aspect of your new Microscope and 2MP Imager. Eight prepared slides are included to get you started. You can also explore the exciting microscopic world by creating your own specimen slides using the included blank slides and tools.

Note: This product is designed and intended for use by those 14 years of age and older.

SPECIFICATIONS

STAGE	Plain Stage with Clips – 90 mm x 97 mm (3.5" x 3.8")
WF 10X AND WF 16X EYEPIECES	Glass Optics
IMAGER	2MP Sensor
SOFTWARE	Digital Microscope Suite - Windows OS: XP, Vista, 7 and 8, Mac OS: 10.4.9 and higher <i>NOTE: No software to install for Mac. Works with Mac based programs such as PhotoBooth and iChat</i>
FOCUSER	Coarse Focus
OBJECTIVES	All Glass Optics – see magnification chart for powers
ILLUMINATOR – TOP AND BOTTOM LED	Uses 4 AA Alkaline batteries (user supplied)
HEAD	Monocular with 45° Incline and 360° Rotatable
NOSEPIECE	Triple with click stop
CONDENSER	N.A. 0.65
WEIGHT / DIMENSIONS (WITH BATTERIES)	3.3 lbs (1.5 kg), 5.25"x8.5"x11.5" (133x216x292 mm)

STANDARD ACCESSORIES WITH YOUR MICROSCOPE

- WF 10x eyepiece
- Digital imager with 2MP sensor
- Imager adapter
- CD ROM with software
- 4x, 10x, 40x (spring loaded) Objective lenses
- Top illuminator – LED with adjustable brightness
- Bottom illuminator – LED with adjustable brightness
- Needle probe and eyedropper
- WF 16x eyepiece
- Filter wheel/diaphragm – 6 positions
- 8 prepared slides and 4 blank slides
- 4 cover slips
- Tweezers

MAGNIFICATION TABLE

Use the following table to determine magnification using your microscope's different eyepiece/objective lens combinations.

NOTE: When using the 2MP Imager, the Imager will act as a 10x eyepiece.

OBJECTIVE LENS	4X	10X	40X
WF 10x Eyepiece	40x	100x	400x
WF 16x Eyepiece	64x	160x	640x

SETTING UP MICRO360

1. Remove styrofoam container from carton.
2. Carefully remove microscope and accessories from container and set on a table, desk or other flat surface.
3. Remove bag covering the microscope.
4. Remove cap from eyepiece tube.
5. Install four AA batteries (user supplied) under the microscope's base (see images below).

6. Insert the WF 10x eyepiece in the eyepiece tube.

MICRO360 OPERATION

Before viewing specimens, please read these sections thoroughly regarding focusing, changing power (magnification), using the stage and adjusting illumination.

ROTATING THE HEAD

Your microscope's head can be rotated 360°. You can view from any position by moving the head (while holding the arm) to a desired location.

VIEWING A SPECIMEN

Carefully place a specimen slide under the stage clips and center the specimen. Images will be observed upside down and reversed right to left, with practice centering objects will become easier to accomplish. Prepared slides, blank slides and tools are included with your microscope to get you started (see images below).

You are now ready to focus and view a specimen. Be careful not to damage a specimen slide or valuable object. When using higher powers while focusing, make sure the objective lens does not hit the slide or object being viewed.

FOCUSING & CHANGING POWER (MAGNIFICATION)

Place specimen slide (or object) directly under the objective lens and use the focus knob to focus on the specimen. For smaller objects, place on a flat material for easier viewing.

1. Always start with lowest power (4x objective lens and WF 10x eyepiece). This objective lens and eyepiece combination will yield 40x – the lowest but most useful power for viewing large or small solid objects.
2. For higher powers, rotate the nosepiece to change the objective lens to 10x or 40x. Hold the microscope above the nosepiece with one hand and rotate the nosepiece with the other hand until it clicks in position. Be cautious not to let the objective lens touch the specimen slide or object when changing to higher powers.

Note: You should first turn the focus knob to lower the stage.

3. You can replace the WF 10x eyepiece with the WF 16x eyepiece to obtain three additional magnifications, including the highest power possible (640x).
4. At higher powers, your views will be greatly magnified but somewhat darker. The most enjoyable views are seen at lower powers which have a wider field of view and brighter illumination.

ILLUMINATION

To achieve the sharpest and best views, proper illumination (lighting) must be used:

1. To power the illuminator(s) on, turn the brightness control switch (located on both sides of base).
2. The top illuminator is used for viewing at low power (4x –objective lens) as higher powers (10x and 40x objective lenses) will block some or all the light. If you need to use high power to observe solid objects, use a bright secondary light (desk lamp, etc.) for direct illumination.
3. The bottom illuminator is used mainly for viewing specimen slides where light shines up through a hole in the middle of the stage through the slide.
4. Having both illuminators on at the same time will provide enough light for thick and irregular specimens.

ADJUSTING THE LIGHTING

Specimens of different sizes, thickness and color variations will require different levels of illumination. You can adjust illumination by turning the brightness control switches as needed.

When viewing a specimen that is not transparent or dark in color, you may need to increase the amount of light to resolve certain features or details. Increase illumination by turning the brightness control switches all the way to their highest setting.

Optimum lighting is found by making adjustments and experimenting. Each specimen may require slightly different illumination as well as same specimens viewed under different magnifications.

When lighting is no longer needed, turn illuminators off to preserve battery life.

USING FILTERS AND DIAPHRAGM

Before using Micro360+, check and make sure no filters are in the optical path. Viewing is normally conducted without colored filters. However, to bring out increased levels of detail, experiment by using different backlighting filter colors for very bright transparent specimens. To change lighting color, rotate the filter wheel to the desired color – Red (R), Green (G), Blue (B). Each color filter is centered once a faint click stop is heard or felt. You may need to adjust the focus knob slightly for best results.

DIAPHRAGM – Within the filter wheel are three diameter openings – (1) 1 mm, (3) 3 mm and (6) 6 mm, which limit the amount of light passing through to the specimen. These openings are part of the diaphragm which allows you to change diameter sizes to maximize contrast, brightness, etc. The default setting is "6".

USING THE 2MP IMAGER

1. Carefully remove the Imager, Adapter, and CD-ROM (Digital Microscope Suite) from the box and set them on a table, desk, or other flat surface.

NOTE: *The USB cable is permanently attached to the Imager.*

2. Remove the plastic bag covering the Imager.

3. Attach the Adapter to the Imager by sliding the Imager into the adapter and tightening the three thumbscrews.

NOTE: *Make sure to first back the thumbscrews out enough, so they do not protrude through inside of the barrel and interfere with the Imager when you insert it.*

NOTE: *The position of the Adapter on the Imager is not critical but the top of the Adapter (thumbscrew end) should be about 40 mm (1.6") from the sensor end for optimal performance on most microscopes.*

4. Remove the eyepiece from the Micro 360+ microscope

5. Put the Imager into the eyepiece barrel. It should slide in easily and be held in place by gravity.

NOTE: *When moving your microscope with the Imager in the eyepiece tube, do not turn the microscope sideways or upside down or the Imager will fall out.*

INSTALL SOFTWARE

MAC:

1. Connect the Imager to your Mac via the USB cable.

2. The Imager is "plug and Play" and the driver should automatically be recognized by your system.

3. Open one of the imaging software programs that come with your Mac, such as Photo Booth or iChat.

4. Your Imager should be recognized by the software and the video stream should be visible.

5. Now adjust the focus with the Micro360+ microscope.

6. Use the software to take images and videos.

NOTE: *Make sure your system is 10.4.9 or higher when using a Mac and at least Windows XP for Windows.*

WINDOWS:

1. Install the CD software on the CD ROM.

2. Connect the Imager via USB.

3. Open the Digital Microscope Suite software.

4. The Imager should be recognized by the software and the video stream should be visible.

5. Use the software to take images and video.

NOTE: *The approximate magnification powers using the imager, are based on the objective lens you are using and the Imager. The Imager magnifies the image approximately 15x. For example, if the objective lens you are using is 4x, then to calculate the total magnification multiply (4x) x (15x)= 60x or 60 times magnification.*

Of course, with digital microscopy, the final magnification also takes into consideration the size of the display that you are viewing on. The bigger the screen, the larger the magnification.

FOCUSING, VIEWING AND CAPTURING IMAGES

1. You will focus using the components of the Micro360+ microscope, in the same way that you would with a traditional eyepiece

2. Viewing is now done on the computer screen

3. To capture an image, focus and then using your computer cursor, select: ACTION- TAKE SNAPSHOT

4.The software is very intuitive and here is a breakdown if each menu item:

- FILE

- New Album: Creates a new album to store images and video
- Print: Allows you to print a selected image in your album
- Email: Allows you to Email a selected image in your album
- File Info: Creates a new album to store images and video
- Settings: Here you can adjust the following:
 - Image/Video Format
 - Image/ Video Resolution
 - Number of photos to take in a photo burst
 - Select your video device (if using multiple devices)
 - Choose the "Media Folder" location on your hard drive for storing images
- Exit: Closes the software

- EDIT

- Copy: Copy a selected image or video file or folder
- Cut: Cut a selected image or video file or folder
- Paste: Paste an image
- Delete: Deletes a selected image or video file or folder

- ACTION

- Take Snapshot- F7: Takes a still image of what is being viewed in the main window
- Record Video- F8: Starts to record a video of what is being viewed in the main window

NOTE: you can also start and stop taking video by pressing the large blue and red button under the main preview window

- Take Multiple Snapshots-F9: Takes multiple still images (based on the number entered in the Settings/Photo Burst window- Default is 5)
- Full Screen Mode-F12: Increases viewing window of preview to full screen. press ESC button to exit full screen

- HELP

- About Digital Microscope Suite 2.0: Gives you general information about the software

- Never disassemble or clean internal optical surfaces. Repair work should be performed by qualified factory technicians or other authorized repair facilities.
- Use care when handling glass specimen slides, as edges can be sharp.

WARRANTY

Your microscope has a two year limited warranty. Please see Celestron's website for detailed information on all Celestron microscopes at www.celestron.com.

2835 Columbia St.
Torrance, California 90503 U.S.A.
©2013 Celestron
All rights reserved. / Printed in China 05-13

This product is designed and intended for use by those 14 years of age and older.

FCC Note: This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

CARE, MAINTENANCE AND WARRANTY

Your Celestron Micro360+ microscope and 2MP Imager are a precision optical instruments and should be treated with care at all times. Follow these care and maintenance suggestions and your microscope will need very little maintenance throughout its lifetime.

- When you are finished using your microscope, remove any specimens left on the stage.
- Turn off top and bottom illuminators when you are finished using your microscope.
- Remove batteries if you will not use your microscope for an extended period of time.
- Store the microscope and imager in a dry and clean place.
- Be very careful if using your microscope and imager in direct sunlight to prevent damage to the microscope and imager or your eyes.
- When moving your microscope, carry it by the "arm".
- Clean outside surfaces with a moist cloth.
- Never clean optical surfaces with cloth or paper towels as they can scratch optical surfaces easily.
- Remove dust with a camel's hair brush or use an air blower.
- To clean fingerprints off optical surfaces, use a lens cleaning agent and lens tissue available at most photo outlets. While cleaning, do not rub in circles as this may cause sleeks and scratches to occur.

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