
Intell-Lab[™] Precision Toploading Balances

PR Series



User Operation Manual

June, 2009

INDEX


1 INSTALLATION INSTRUCTIONS	1
2 WEIGHING PAN ASSEMBLY	2
3 KEYBOARD AND DISPLAY	3
4 INPUT AND OUTPUT CONNECTORS.....	4
4.1 REAR AND BOTTOM PART OF BALANCE, PR MODEL	4
5 STANDARD WEIGHING	5
5.1 STAND BY.....	5
5.2 SIMPLE WEIGHING	5
6 CALIBRATION	6
6.1 EXTERNAL CALIBRATION BALANCES	6
6.1.1 <i>External calibration</i>	6
6.2 BALANCES WITH INTERNAL CALIBRATION	8
6.2.1 <i>Internal calibration</i>	8
6.2.2 <i>Autocalibration (A UT-CAL)</i>	8
6.2.3 <i>Internal calibration (I-CAL)</i>	9
6.2.4 <i>External calibration (E-CAL)</i>	9
6.2.5 <i>Technical calibration (TEC-CAL)</i>	9
7 TARE FUNCTION.....	11
7.1 OPTIONAL ALPHANUMERIC EXTERNAL KEYBOARD (CODE T20 1)	12
7.2 MANUAL TARE FUNCTION	13
8 WEIGHT UNITS	14
9 PC COMMUNICATION SETTING.....	16
10 PRINTER SERIAL COMMUNICATION SELECTION	17
11 TRANSMISSION SPEED SELECTION.....	18
12 AUTOZERO FUNCTION.....	19
13 FILTERS SELECTION	20
14 STABILITY FUNCTION	21
15 BACKLIGHT SETUP.....	22
16 AUTO POWER-OFF FUNCTION	23
17 PIECECOUNTING FUNCTION	24
17.1 VISUALIZATION OF TOTAL AND UNIT WEIGHT OF PIECES.....	25
17.2 AUTOMATIC UPDATE OF UNIT WEIGHT	26
17.3 MANUAL INSERTION OF THE UNIT AVERAGE WEIGHT.....	27
18 V-RANGE FUNCTION	28
19 DENSITY DETERMINATION OF A SOLID OR A LIQUID	29
19.1.1 <i>DENSITY DETERMINATION OF A SOLID</i>	29
19.1.2 <i>DENSITY DETERMINATION OF A LIQUID</i>	30
20 MAXIMUM LOAD DETERMINATION FUNCTION	32
21 RS232 INTERFACE FEATURES	33

22 CONNECTORS PLACES (REAR).....	37
23 WORKING WITH BATTERY.....	38
24 ERROR CODES	39
25 QUICK GUIDE TO BALANCE PARAMETERS SETUP.....	40
26 QUICK GUIDE TO THE USE OF THE BALANCE PROGRAMS.....	41
27 BALANCE TECHNICAL CHARACHTERISTICS	42


1 Installation instructions

- ☐ **Remove** balance and accessories from the cartoon.
- ☐ **Do not install** the balance in a place with air flows, heavy thermic changes and vibrations.
- ☐ **Place** the weighing pan and the support pan on the balance (pag. 2-3).
- ☐ **Level** the balance using the level bubble and levelling feet located underneath the case
- ☐ **Connect** power supply to connector 2 located on the rear panel of the unit (page 5).
- ☐ **Insert** power cord into AC outlet, **which shall be easily accessible**; after few seconds the balance will automatically switch on.
- ☐ **Wait 30 minutes from switch on** and then calibrate the balance using the mass enclosed in the package, following the instructions (page 7-9).
- ☐ **Calibrate** the balance again every time it is removed from previous place.
- ☐ **Check** balance calibration periodically.
- ☐ **We recommend** not to drop heavy objects on balance pan, in order to avoid damages.
- ☐ **Assistance service** must be effected by specialized staff and the spare parts used must be original. Therefore, it is necessary to apply to the reseller who sold the equipment.
- ☐ **Keep package** in case the balance should be sent back to the factory for service.


2 Weighing pan assembly

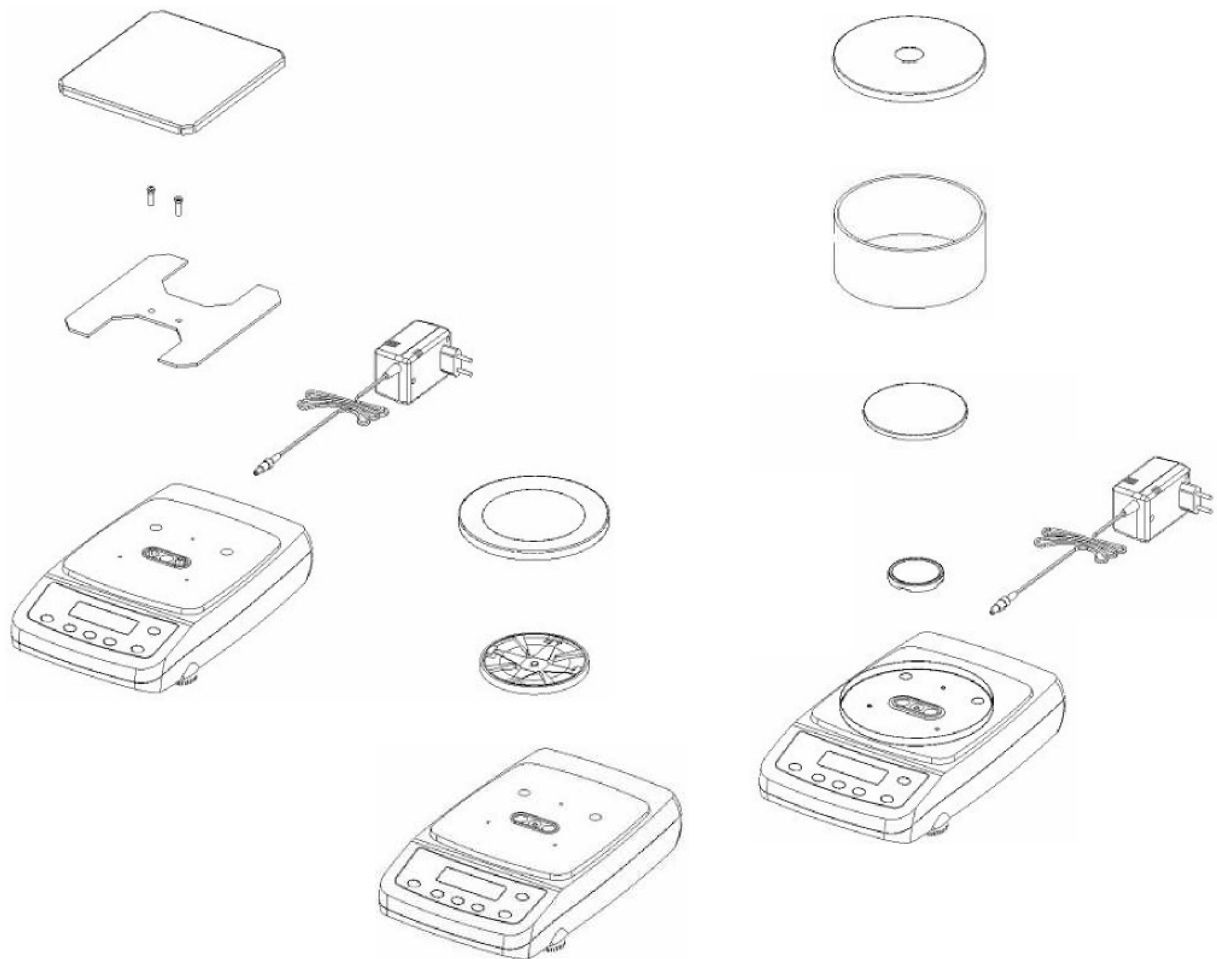


3 Keyboard and display

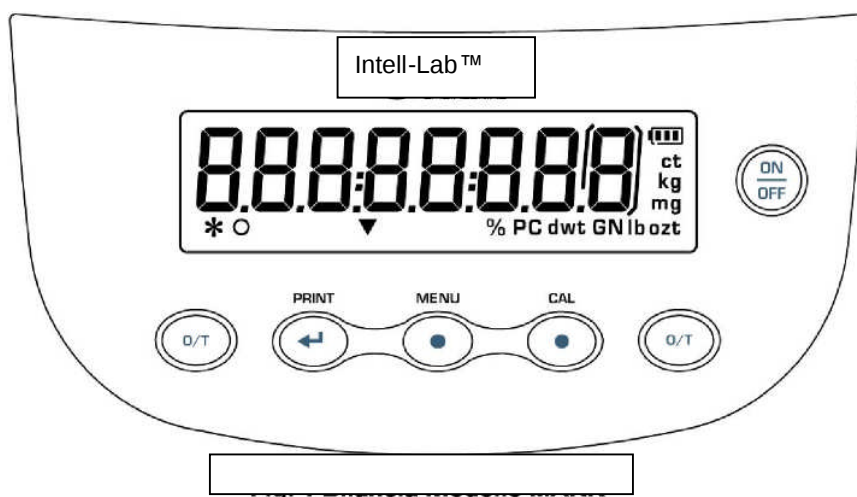


Standby (OFF) or power on (ON).

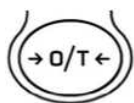


TARE or zero button.

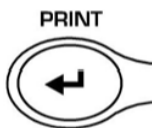

Selection CONFIRM or SEND data to printer.

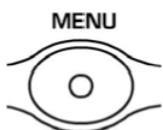


Balance setup MENU button.

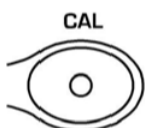

Balance CALIBRATION button.

4 Input and output connectors

4.1 Rear and bottom part of balance, L model

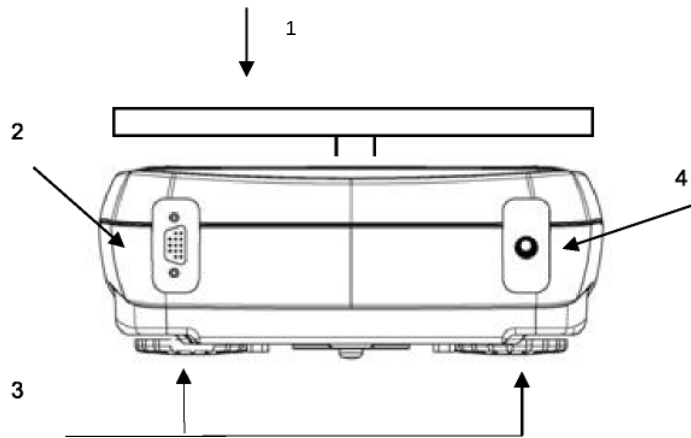


1. Weighing pan

2. **CONNECTOR 1**

9 pin RS232 Interface
Output for PC/keyboard or
printer

3. Levelling feet

5. **CONNECTOR 2** power
supply connector

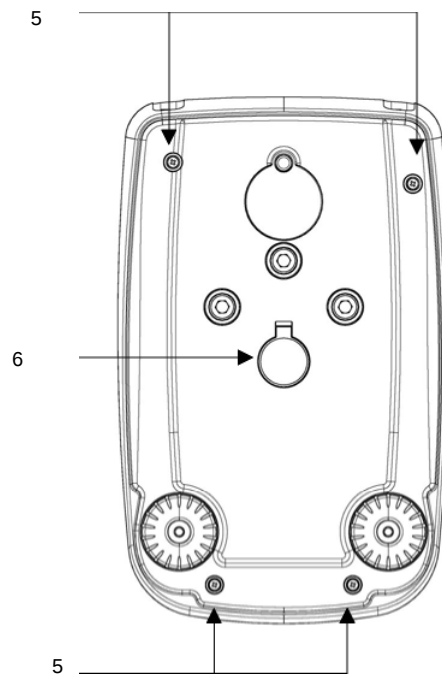


5. Screws for balance closing

6. Hook for weighing under
balance pan. First remove
hook cap.

5 Standard weighing

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
After having connected the balance to AC outlet, it will perform an internal circuits test, after that the balance will set itself in stand-by mode.



5.1 Stand By


From “**STAND BY**” mode:

- ☐ Press **ON/OFF** button to bring balance to work conditions.
- ☐ Press again **ON/OFF** button to return to “**STAND BY**” condition.

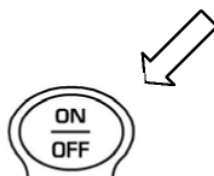

5.2 Simple weighing

Load the sample to weigh on the pan and read the value on display as soon as the stability symbol (*star*) appears.

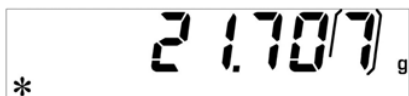

6 Calibration

Electronic balances take mass measurements making use of gravity (g). Difference of latitude in geographic areas and altitude will vary gravity acceleration value (g).


Therefore, for accurate measurements, the balance must be adjusted to the local environment. This adjustment is accomplished by calibration function.

6.1 External calibration balances

6.1.1 External calibration

Calibration is accomplished by pressing **CAL** button.

1. Press **CAL** button when pan is empty, dashes are displayed on the display..

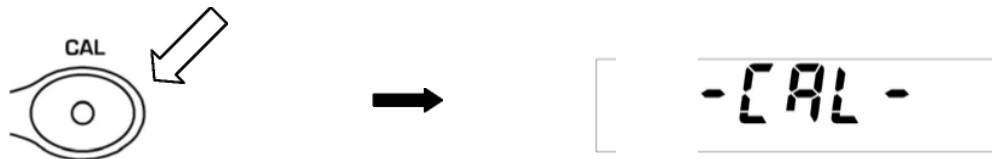

2. When calibration weight value starts to flash, load the weight on the pan.




The display will stop flashing, indicating calibration weight value. Wait for the acoustic signal to confirm calibration.

3. Unload calibration weight from the pan. The balance is ready for weighing operations.

NOTE: if there is an interference during calibration process, an error message will be displayed.


Moreover, it is possible to calibrate the balance with a calibration weight higher than the one set by default:

1. Press and keep **CAL** button pressed with empty pan until the acoustic alarm stops, then release the button. On display it will be visualized the string "-CAL-", followed by flashing string "**LOAD**".

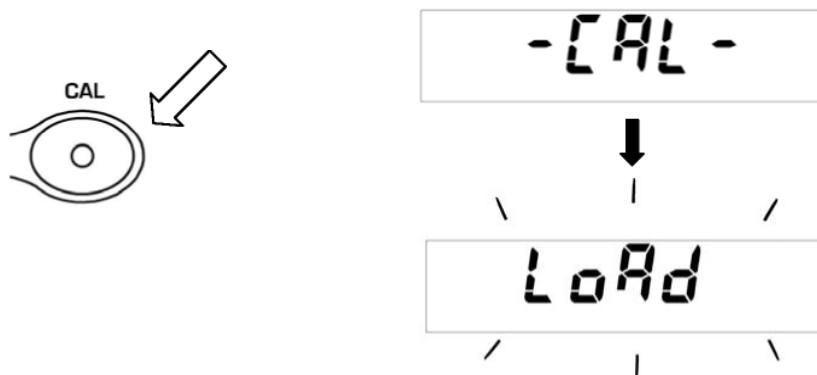


2. Load on the pan a weight equal or higher than default calibration weight; the balance will recognize it as valid weight if equal or higher than calibration weight as long as it is a whole number in comparison with the most meaningful digit of calibration weight . e.g.: *if calibration weight is 200g, it will be possible to calibrate the balance with values from 200g, 300g, 400g up to the highest limit of balance weighing range.*

The message "**LOAD**" on display will stop flashing. Once calibration has been effected, the value of calibrated weight will be displayed.

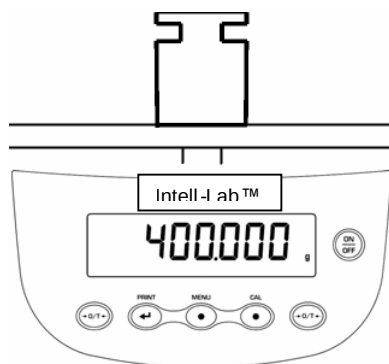

3. Unload calibration weight. The balance is ready for weighing operations.

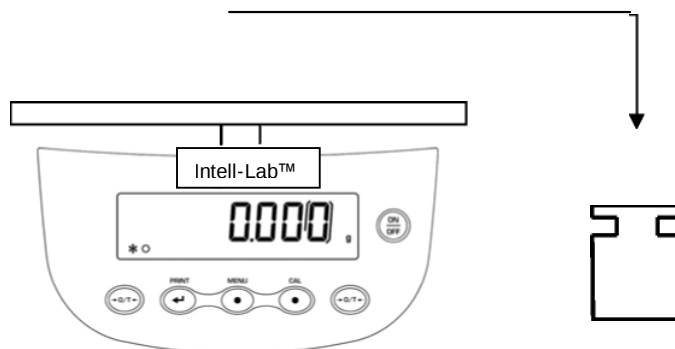

NOTE: if there is an interference during calibration process, an error message will be displayed.

6.2 Balances with internal calibration


6.2.1 Internal calibration

In these balance models there are 4 calibration modes:

From display zero condition, press and keep pressed the **MENU** button until the acoustic alarm is over, then release the button. The message “**units**” will be visualized on display, press then **MENU** button until you visualize “**Calib**” on display. Press ENTER to confirm.

1. Select the calibration mode you wish by pressing **MENU** button in sequence:

- ~ **AUT-CAL**: autocalibration
- ~ **I-CAL**: internal calibration
- ~ **E-CAL**: external calibration
- ~ **TEC-CAL**: technical calibration

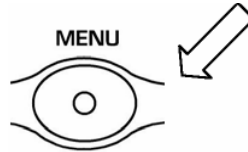



2. Press **PRINT** button to confirm “**AUT-CAL**”, “**I-CAL**”, “**E-CAL**”. To confirm “**TEC-CAL**” keep pressed the **PRINT** button until the acoustic alarm is over.
3. After selection, the balance returns to calibration menu. Press and keep pressed **MENU** button until the acoustic alarm is over, then release the button. Balance is again ready for weighing operations.

6.2.2 Autocalibration (AUT-CAL)

The balance autocalibrates through the internal reference calibration mass, after the microprocessor has checked that no other weighing operations are being effected. In this mode it is furtherly possible to effect calibration through reference internal mass by pressing **CAL** button at any time, being sure first that there's no weight load on the pan.

1. Press **CAL** key at empty pan.

The message “**CAL**” will be displayed and balance calibration will be effected automatically.

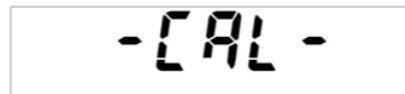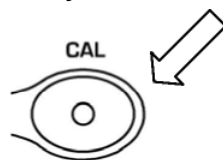

2. At the end of calibration, balace returns to normal weighing conditions.



If calibration is not finished due to vibrations or air flows, then the message “**CAL bUt**” is displayed. Press again **CAL** button. If the problem repeats then select external calibration and contact the product supplier.

6.2.3 Internal calibration (I-CAL)

The balance calibrates itself with internal reference mass **ONLY** on user command by pressing **CAL** button.

Before effecting internal calibration, be sure that there's no weight loaded on the balance pan.

6.2.4 External calibration (E-CAL)

The balance is calibrated through a external reference mass (**NOT** provided with models with internal calibration).

(Follow procedures described at paragraph 6.1.1)


6.2.5 Technical calibration (TEC-CAL)

This function allows to store the value of internal reference mass whenever checking or assistance actions require it.


1. After having selected the **TEC-CAL** calibration mode, press **CAL** button at empty pan. It will be displayed "CAL".




2. When the value of calibration weight start flashing on display, load the weight on to the balance pan.

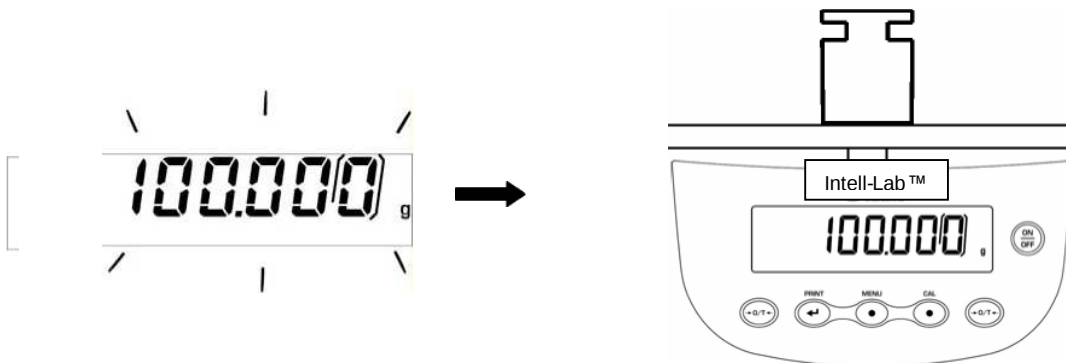

3. Wait the acoustic alarm and that the displayed calibrated weight value stops flashing, then unload the weight from balance pan.

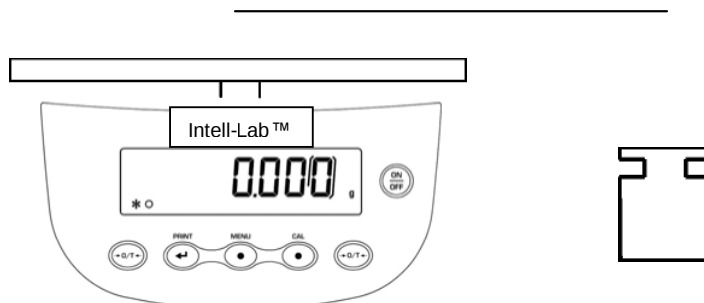

4. When string “0.000” is displayed continuously, then press and keep pressed the **PRINT** button. This starts the internal weight value automatic acquisition and store. During the acquisition cycle, the display will show “tEC-MEM”.

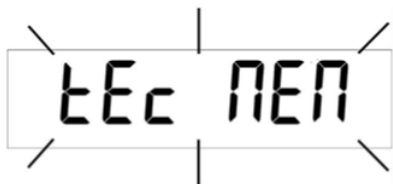

5. Once stored the value of internal calibration weight, balance returns to normal weighing conditions.
6. Return to calibration menu as described at paragraph 6.2 and set the desired calibration mode: internal, automatic or external.

ATTENTION : this procedure must be effected only using E2-class reference masses.

7 Tare function

1. Load the container on the pan. The display will show the weight.




2. Press **O/T** button. “**O-t**” string will be displayed.



3. After reaching stability, the value “**0.000**” will be displayed. If the stability is not reached (due to air flows or vibrations or other disturbs) the dashes will remain displayed.



4. Load the objects to weigh in the container. Read net weight value on display.

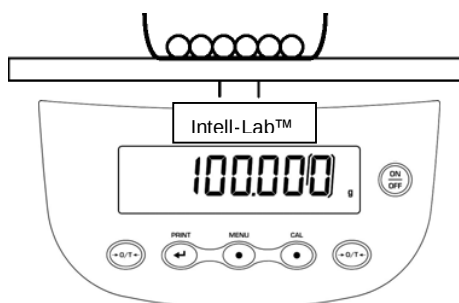

7.1 Optional alphanumeric external keyboard (code T201)


If you have the optional external alphanumeric keyboard, then it is possible to perform tare also by pressing the **TARE/DELETE** button of this keyboard, in the same way as previously described.

UP SELECT	TARE DELETE	TARE HIV. M		HELP STAT.	
DOWN FORMUL	CLEAR	ESCAPE		INSER	
A	B	C	D	E	
F	G	H	I	J	
K	L	M	N	O	
P	Q	R	S	T	
U	V	7	8	9	
W	X	4	5	6	
Y	Z	1	2	3	
/	-	.	0	SP	

7.2 Manual tare function

If you have the optional external alphanumeric keyboard (T201), it is possible to insert manually a known value of tare.

1. Press **TARE MAN** button on the alphanumeric keyboard. An arrow will appear on the balance display along with the previous value of manual tare, if one has been inserted before.

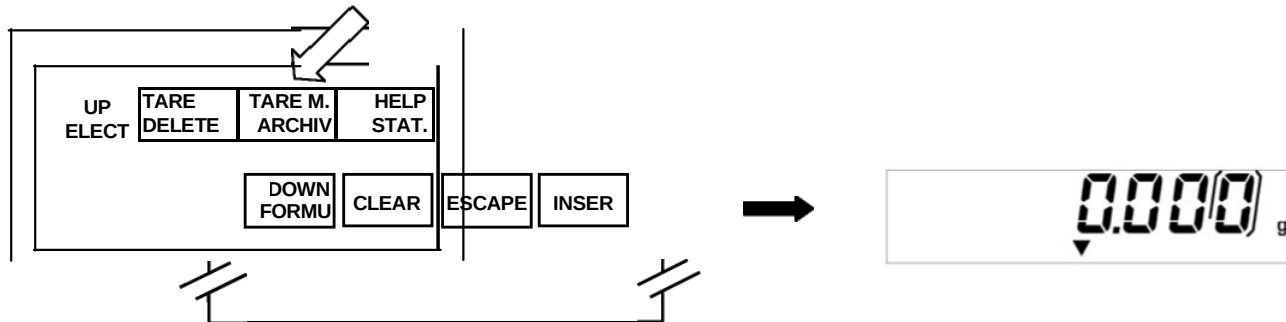

2. Press **CLEAR** to set at zero the previous value (if any) then insert the new value using the alphanumeric buttons placed on the bottom part of alphanumeric keyboard.



3. Press **INSERT** to confirm.
4. Press **ESCAPE** button to escape from tare condition.

To set to zero the value of the manual tare inserted by means of alphanumeric keyboard, effect a normal operation of tare by pressing **O/T** button of the balance or with **TARE/DELETE** button placed on the optional alphanumeric keyboard

NOTE: manual tare function is **ONLY** available for models equipped with alphanumeric keyboard.

8 Weight units

It is possible to select the weight unit that balance will use to display weight.


1. From display zero condition press and keep pressed the **MENU** button until the acoustic alarm is over, then release the button. The message “unitS” will be displayed then press **PRINT** to confirm..
2. It will



displayed “GRAM” unit. Pressing now the **MENU** or **CAL** button, it will be possible to scroll forward or backward the weight units menu.



3. Premere il tasto **PRINT** per confermare o **MENU** per passare ad un'altra unità di misura.

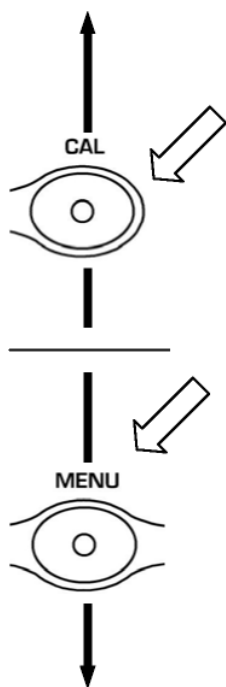

SYMBOL	WEIGHT UNIT	CONVERSION FACTOR 1g =
GRAM	GRAMS	1.
CARAT	CARATS	5.
OUNCE	OUNCE	0.035273962
POUND	POUNDS	0.0022046226
PENN.	PENNYWEIGHTS	0.643014931
ONCETR.	OUNCE TROY	0.032150747
GRANO	GRAIN	15.43235835
Tael HON	HONG KONG Tael	0.02671 725
* Tael SGP	SYNGAPORE Tael	0.02646063
* Tael ROC	R.O.C. Tael	0.02666666
* MOMME	MOMME	0.2667
* PER 10	x 10	10.
* PER 100	x 100	100.

* Not available for all models

4. After having selected the desired weight unit, press the **MENU** button to go to next parameter or **CAL** button to return to previous.
5. To escape from parameters setup menu, press the **MENU** button until the acoustic alarm is over, then release the button. The balance returns to normal weighing conditions.

9 PC communication setting

Connect the balance to PC with proper cable (pag. 35)

1. From zero display condition press and keep pressed the **MENU** button until the acoustic alarm is over, then release the button. It will be displayed the string "**units**", then press the **MENU** button until it is displayed "**PC-PRTR**" message then confirm by pressing **PRINT** button.

2. Press button

"PC


to

print or

to



print to

command, then press the **PRINT** button to confirm the choice.

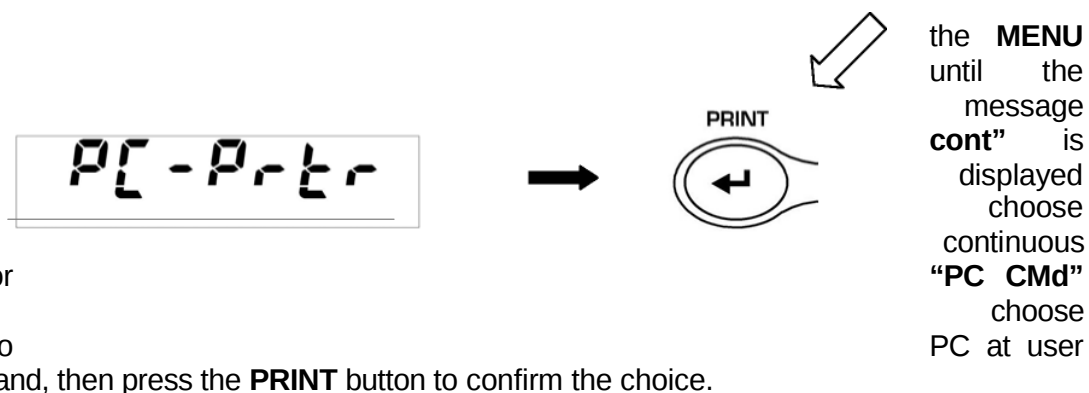

PC Cont

PC Cmd

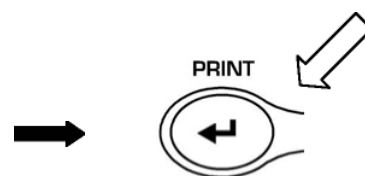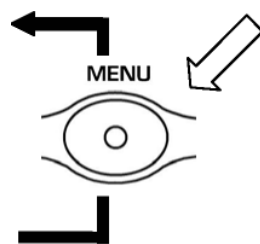

3. After having selected the desired transmission mode, press the **MENU** button to go to next parameter or the **CAL** button to go to previous.
4. To escape from parameters setup menu, press the **MENU** button until the acoustic alarm is over, then release the button. The balance returns to normal weighing conditions.
5. The balance returns to normal weighing conditions sending the data continuously or at user command.





NOTE: select speed transmission (pag. 19)

10 Printer serial communication selection

Connect the balance to the printer using the proper cable (pag. 35)




1. From zero condition on display, press and keep pressed the **MENU** button until the acoustic alarm is over, then release the button. The message “**unitS**” will be displayed, then press **MENU** until you read the message “**PC-PRTR**” on display and confirm pressing **PRINT** button

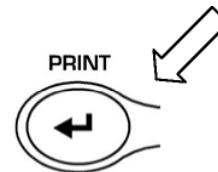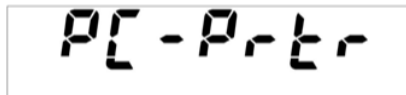

2. To select the data print mode, press

MENU until the message “**PRINT**” is displayed.

3. Press **PRINT** to confirm.

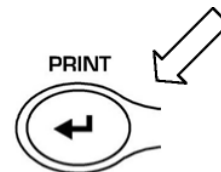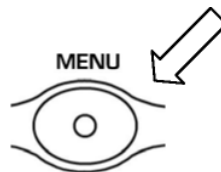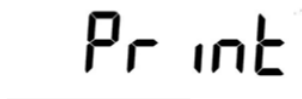

4. After having selected the data print mode, press **MENU** button to go to next parameter or **CAL** button to go to previous one.
5. To escape from parameters setup menu, press **MENU** button until the acoustic alarm is over, then release the button. The balance returns to normal weighing conditions, ready to transmit data each time the **PRINT** button is pressed.




NOTE: select speed transmission (pag. 18)

11 Transmission speed selection

1. From zero condition on display, press and keep pressed the **MENU** button until the acoustic alarm is over, then release the button. The message “unitS” is displayed, then press **MENU** button until the message “BAUD RT” is displayed and confirm by pressing the **PRINT** button.

baud rt

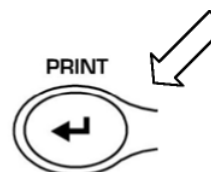

2. Select serial data

transmission speed (1200-2400-4800-9600 baud). Pressing **MENU** or **CAL** buttons it will be possible to scroll forward or backward through the different transmission speeds, then confirm your choice by pressing **PRINT** button.

br 1200






br 2400



br 4800



br 9600

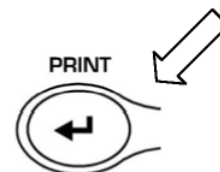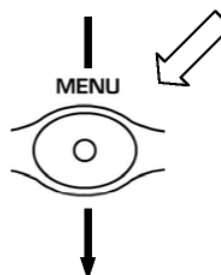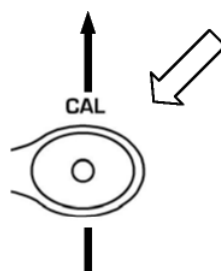

3. After having selected the transmission speed you wish, press the **MENU** button to go to next parameter or **CAL** button to go to previous one.
4. To escape from parameters setup menu, press the **MENU** button until the acoustic alarm is over, then release the button.
5. The balance will return to standard weighing conditions.

*O 0.0000 g

12 Autozero function


Autozero is an automatic correction of a possible zero drift.

- ~ **Au0 OFF** = autozero disabled
- ~ **Au0 1** = light autozero
- ~ **Au0 2** = medium autozero
- ~ **Au0 3** = heavy autozero
- ~ **Au0 3E** = heavy autozero over all range

1. From zero condition on display, press and keep pressed the **MENU** button until the acoustic alarm is over, then release the button. The message “**units**” is displayed, then press **MENU** button until the message “**AUTO 0**” is displayed, then press **PRINT** to confirm.

 →

2. Pressing **MENU** or **CAL** button it will be possible to scroll forward or backward through the different autozero levels, select the one you wish and confirm it by pressing the **PRINT** button.

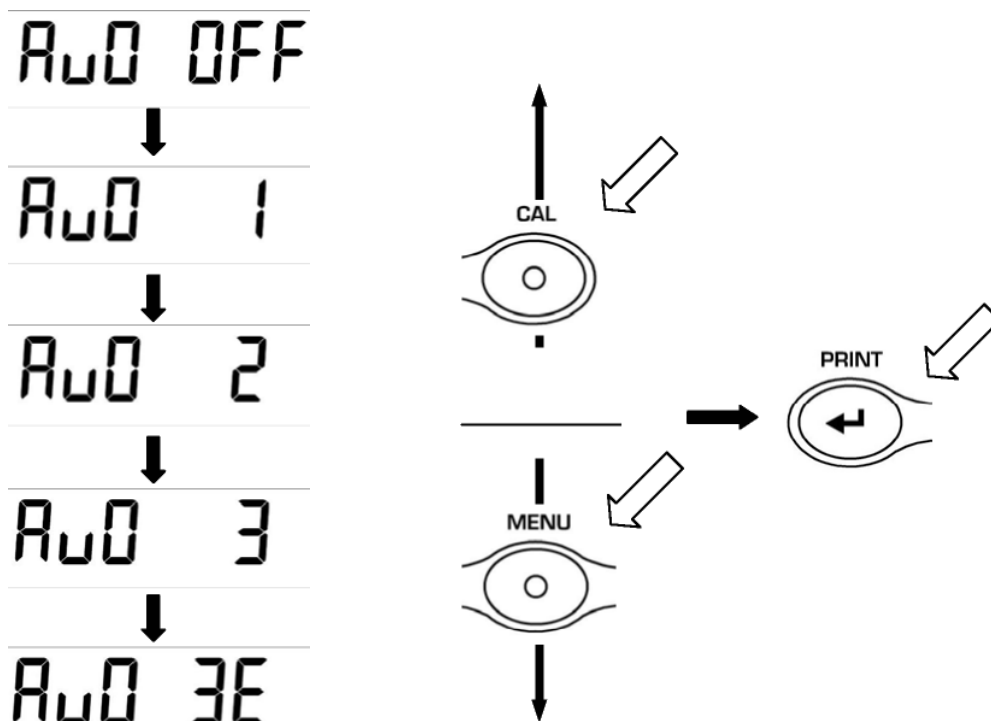

3. After having selected the autozero level you wish, press the **MENU** button to go to next parameter or **CAL** button to go to previous one.
4. To escape from parameters setup menu, press the **MENU** button until the acoustic alarm is over, then release the button.
5. The balance will return to standard weighing conditions.

13 Filters selection


It is possible to adapt the balance to the different environment conditions thanks to the selection of three filters:

- ☐ **FILTER 1:** proportion of ingredients
- ☐ **FILTER 2:** stable conditions
- ☐ **FILTER 3:** unstable conditions

1. From zero condition on display, press and keep pressed the **MENU** button until the acoustic alarm is over, then release the button. The message “**units**” is displayed, then press **MENU** button until the message “**FILTER**” is displayed then confirm it by pressing the **PRINT** button.



2. Pressing **MENU** or **CAL** button it will be possible to scroll forward or backward the different filtering levels, select the one you wish and then confirm it by pressing the **PRINT** button.

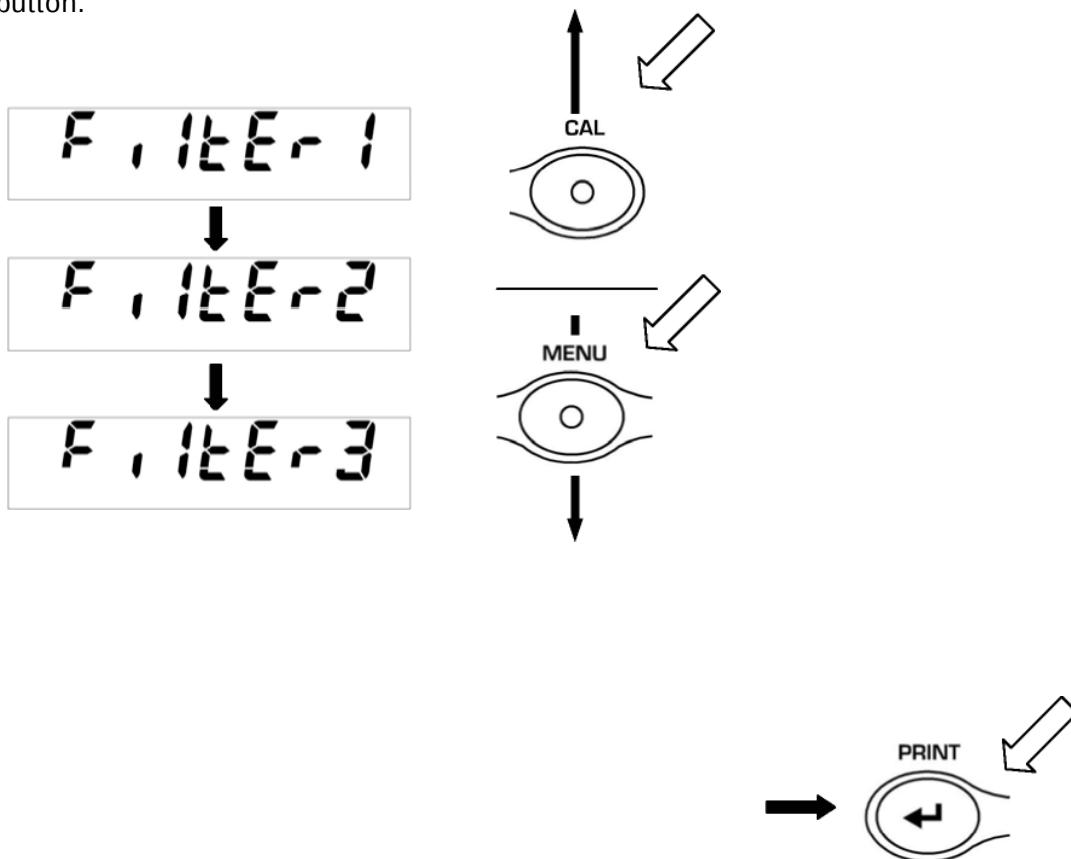

3. After having selected the filtering level you wish, press the **MENU** button to go to next

parameter or **CAL** to go to previous one.

4. To escape from parameters setup menu, press the **MENU** button until the acoustic alarm is over, then release the button.
5. The balance will return to standard weighing conditions.



NOTE: It is suggested to use FILTER 1 when proportion of ingredients must be performed

1. The balance returns to normal weighing conditions.

14 Stability function


The stability symbol is displayed when the weight is stable inside a defined range.

- ☐ **STAB 1** = for stable environments
- ☐ **STAB 2** = for not so stable environments
- ☐ **STAB 3** = for unstable environments


1. From zero condition on display, press and keep pressed the **MENU** button until the acoustic alarm is over, then release the button. The message “**units**” is displayed, then press **MENU** button until the message “**StAbiL**” is displayed, then confirm this by pressing the **PRINT** button.



2. Pressing **MENU** or **CAL** button it will be possible to scroll forward or backward the different stability levels, select the one you wish and then confirm it by pressing the **PRINT** button.

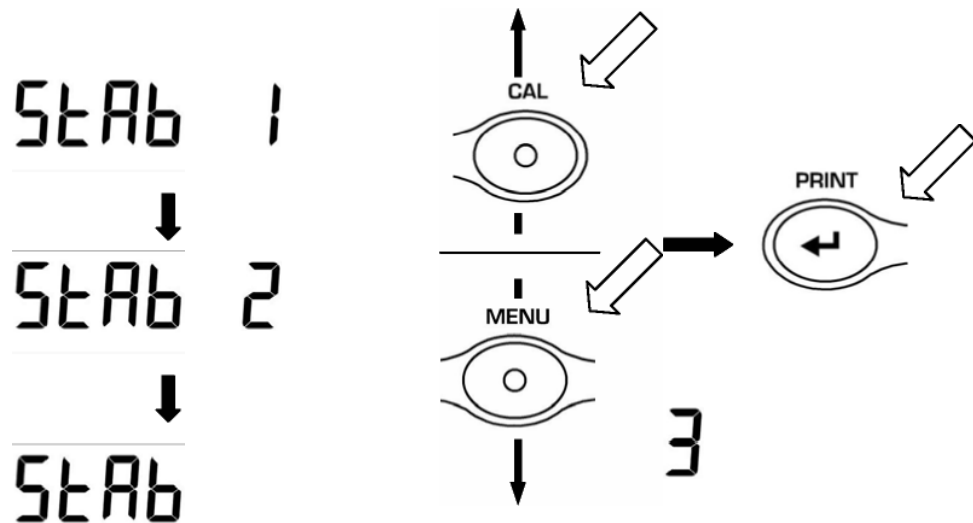

3. After having selected the stability level you wish, press the **MENU** button to go to next parameter or the **CAL** button to go to previous one.
4. To escape from parameters setup menu, press the **MENU** button until the acoustic alarm is over, then release the button.

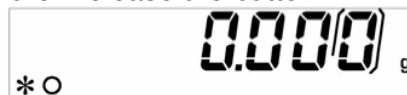


15 Backlight setup

The balance display is equipped with backlight to make the indication more visible also during low light conditions.


There are 3 working modes:

- ☐ **ON** = backlight always switched ON
- ☐ **OFF** = backlight always switched OFF
- ☐ **AUTO** = backlight automatically switched on during weighing operations

1. From zero condition on display, press and keep pressed the **MENU** button until the acoustic alarm is over, then release the button. The message “unitS” is displayed, then press **MENU** button until the message “bLt ”, then press the **PRINT** button to confirm this.

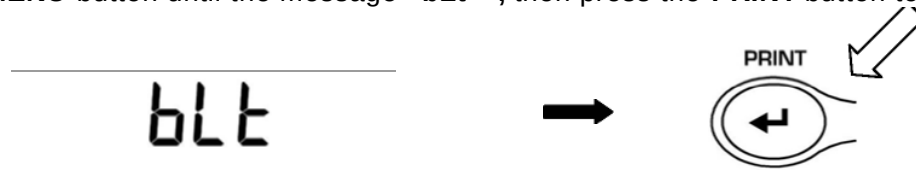

2. Pressing **MENU** or **CAL** button it will be possible to scroll forward or backward the different working modes, select the one you wish and then confirm it by pressing the **PRINT** button

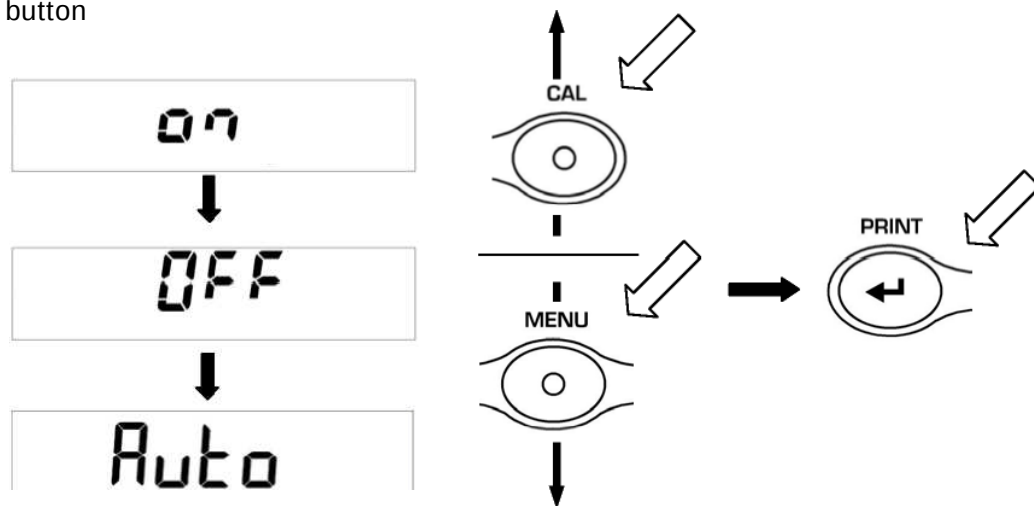

3. After having selected the backlight working mode you wish, press the **MENU** button to go to next parameter or the **CAL** button to go to previous one.
4. To escape from parameters setup menu, press the **MENU** button until the acoustic alarm is over, then release the button.
5. The balance returns to normal weighing conditions.

1. The balance returns to normal weighing conditions.




16 Auto Power-off function

This function allows to activate the automatic power-off of the balance after a defined balance idle-time.

There are 4 auto power-off modes:



- ~ **disab** = Auto power-off disabled
- ~ **2 Min** = Auto power-off after 2 minutes of idle time
- ~ **5 Min** = Auto power-off after 5 minutes of idle time
- ~ **15 Min** = Auto power-off after 15 minutes of idle time

- From zero condition on display, press and keep pressed the **MENU** button until the acoustic alarm is over, then release the button. The message “**units**” is displayed, then press **MENU** button until the message “**time off**” then confirm by pressing the **PRINT** button.

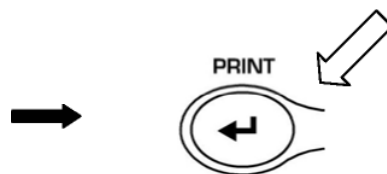

- Pressing **MENU** or **CAL** button it will be possible to scroll forward or backward the different auto power-off modes, select the one you wish and then confirm it by pressing the **PRINT** button

- After having selected the auto power-off mode you wish, press the **MENU** button to go to next parameter or the **CAL** button to go to previous one.

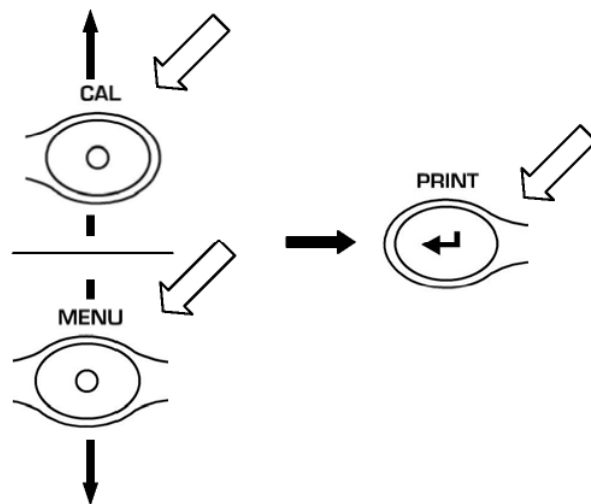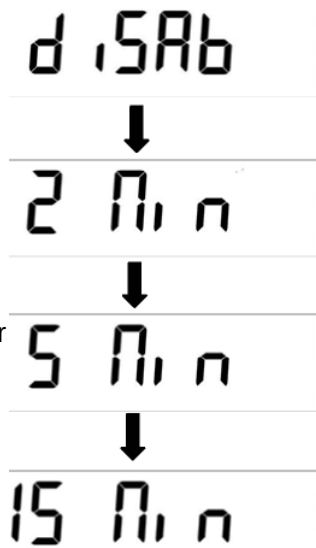


- To escape from parameters setup menu, press the **MENU** button until the acoustic alarm is over, then release the button.



17 Piececounting function

** Not available for all models **


1. From zero condition on display, press **MENU** button until the message “**Count** “ is displayed then press the **PRINT** button to confirm

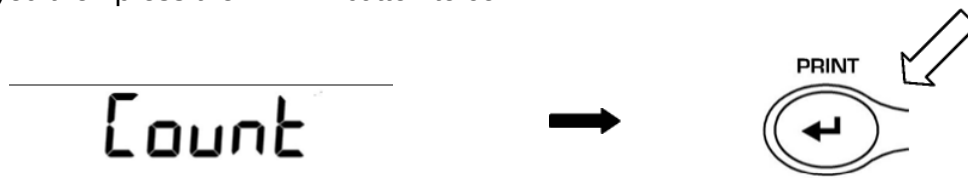

2. Select the number of pieces to put on the pan as sample by pressing **MENU** repeatedly to increase and **CAL** to decrease the number. Choice of the number of pieces (10, 25, 50, 100) depends from the weight of single pieces.
If available, load the empty container first then press **PRINT** to confirm.




Load on the pan the number of pieces displayed on the balance display

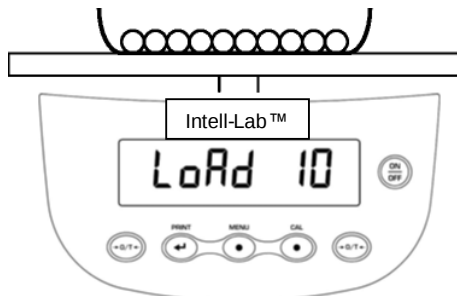


Press again the **PRINT** button and wait the weight to stabilize.

If the number of samples is enough (for example 10 as in figure), then this number will be displayed and it will be possible to go on by loading the pieces to count on the balance pan.

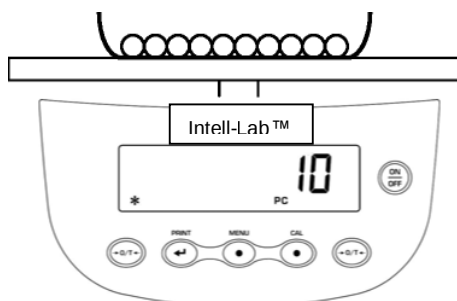


If the weight of the pieces to count is too low in comparison with balance resolution, then an error message will be displayed. In this case it is necessary to use a balance with higher resolution.

If the weight of samples is acceptable but not enough, then the message “**Add**” will be displayed.
Add a quantity of pieces so that to almost double the number read on display, then press the **PRINT** button.

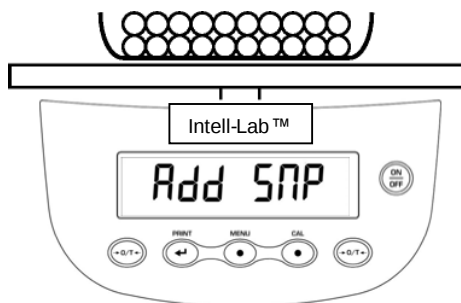

If the number of pieces is still not enough then the message “**Add SMP**” is displayed. Add again a quantity of pieces so that to almost double the number you read on display


As soon as it is reached a sufficient number of pieces, this number is displayed and it will be possible to go on with count by loading the pieces to count on the balance pan.

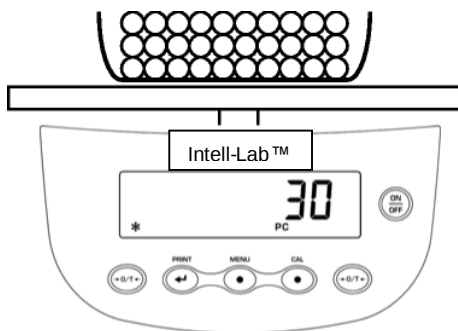

3. To escape from piececounting mode press the **ON/OFF** button and the balance returns to normal weighing conditions.

17.1 Visualization of total and unit weight of pieces

1. Press **MENU** button to pass from pieces number visualization to total weight visualization.

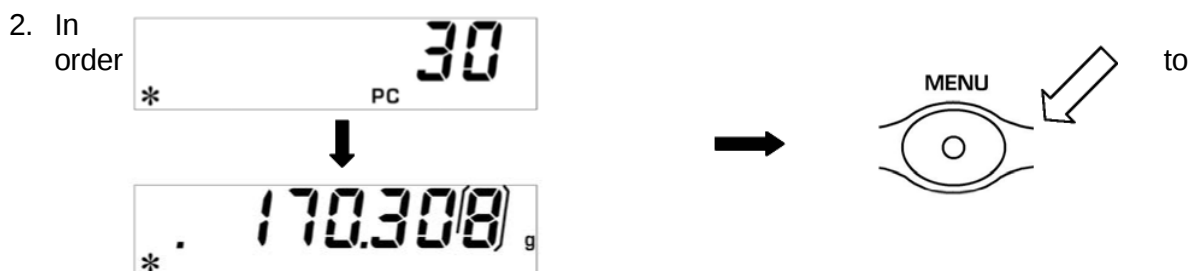


display the weight of single piece from the total number of pieces press and keep pressed the **MENU** button until the acoustic alarm is over.

3. Press again the **MENU** button to display the number of pieces.


17.2 Automatic update of unit weight

After having effected the sampling, it is possible to update the average piece weight as follows:

1. Instead of loading all the pieces to count, load a number of pieces equal to the double of the ones loaded on the pan and wait for the acoustic alarm.

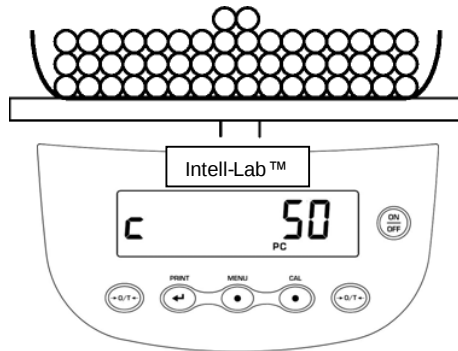

2. It is now possible to repeat this procedure upto a maximum of 255 pieces or to proceed to the normal counting of the pieces.
This procedure allows a more accurate evaluation of the average unit weight and a better precision in the pieces count.

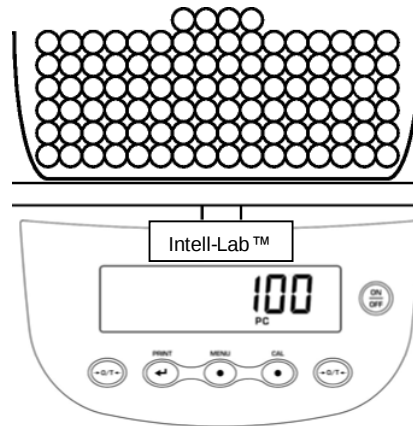

NOTE: automatic update is not active if the sampling has been effected by insertion of the average unit weight from alphanumeric keyboard.


17.3 Manual insertion of the unit average weight

It is possible to use the optional alphanumeric keyboard to insert the weight of the unit sample, if known.

1. From zero condition on display, press the **MENU** button until the message "**Count**", then press the **PRINT** button to confirm.

Select **MANUAL** by pressing repeatedly the **MENU** button.

2. Press button to

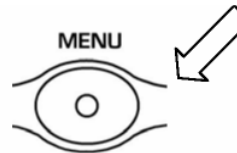

the **PRINT** confirm.

3. Insert

the unit weight of sample in

a

grams using the numerical buttons from 0 to 9 and the decimal point.

In case of mistake, press the **CLEAR** button to start insertion again.

Press the **INSER** button to confirm.

If the inserted weight is 100 times lower than balance resolution, it will be displayed an error message.

To escape without inserting the weight press **ESCAPE** (on alphanumeric keyboard) or **ON/OFF**.


4. If the weight is enough then "**0**" is displayed ; it is now possible to go on to count loading the pieces on the pan.
5. Press **MENU** button to display the total number of pieces, then press again the button to display again the number of pieces.
6. To escape from piececounting function, press **ON/OFF** button.

NOTE: Available **ONLY** for models provided with alphanumeric keyboard (T201).

18 V-Range function

* This (optional) function is available ONLY for double scale models *

1. V-RANGE function is to be considered when it is necessary to effect proportioning of small quantities with accuracy. To activate **V-RANGE** function, after having put on the pan the container if any, press and keep pressed the **TARE** button until the acoustic alarm is over. In this way a tare operation is automatically performed, at the end of which it is possible to operate with the highest resolution range.

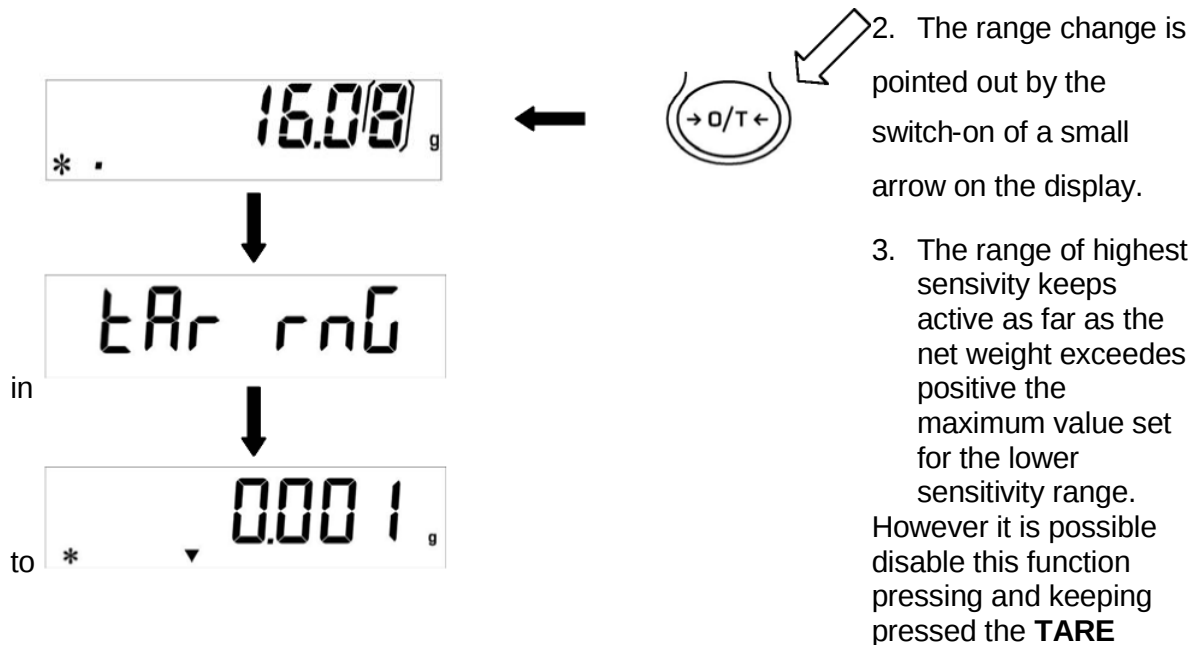

button until the acoustic alarm is over. In both the cases total weight is displayed.

4. In V-Range function the manual tare insertion by (optional) alphanumeric keyboard is not available.

19 Density determination of a solid or a liquid


* The balance must be equipped with the proper optional software (not available for all models) *

The software for density determination allows to calculate the density of a solid or a liquid through the under balance hook or through the optional hydrostatic kit (T002 / T027)

19.1.1 DENSITY DETERMINATION OF A SOLID


1. From zero condition on display, press the **MENU** button until the message “**dEnS**” is displayed, then press the **PRINT** button to confirm.

dEnS

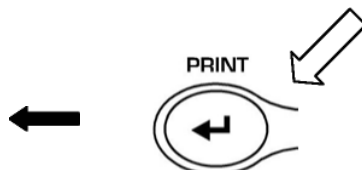


2. Select the function “**d SoLid**” through the MENU button and then confirm with **PRINT** button.

dSoL id

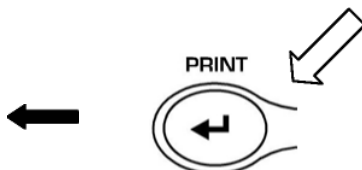

3. The density of the liquid to use will be displayed, the default value is equal to 1.0000 (distilled water at 20°C). It is possible to set a different value if the balance is equipped with the optional alphanumeric keyboard. Once the wished value is set, press **PRINT** button.


dL1.0000

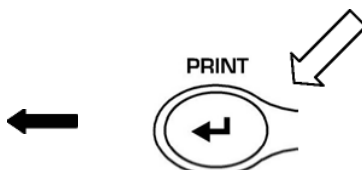

4. Now it will be asked to weigh the solid in air, confirm this with **PRINT** button.

5. If necessary, operation and solid. Press button to value. During acquisition the **WEI AIR** will be flashing.


WEI AIR

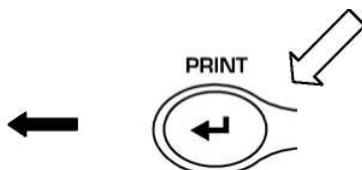

effect a tare load the the **PRINT** acquire the value message

6. Now it is asked to weigh the solid inside the liquid. Effect the tare of the small basket inside liquid, immerse the solid and press **PRINT**. During the acquisition of the value the message **WEI LIq** will be flashing.

WEI LIq

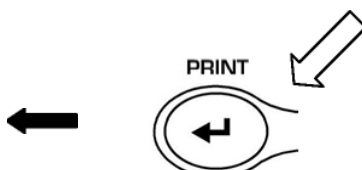

- The result of density of the solid is then displayed. Pressing the **PRINT** button it is possible to print the value of density, if the balance is equipped with a printer.



- If any error occurs, this message will be displayed:




- Press now the **ON/OFF** button to escape from density function or the **MENU** button to perform another measurement.

19.1.2 DENSITY DETERMINATION OF A LIQUID

- From zero condition on display, press the **MENU** button until the message “**dEnS**” is displayed, then press the **PRINT** button to confirm.


dEnS

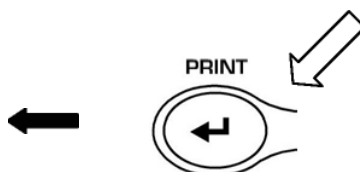

- Using the **MENU** button, select the function “**d Liquid**” and confirm this with **PRINT**

- It will be the value density of float of volume to default

dL 1901d

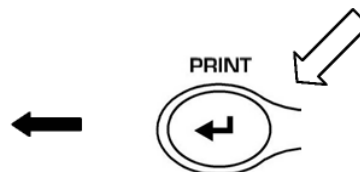


displayed of the the glass known use, the value is

equal to 3.0000. It is possible to set a different value if you have the optional alphanumeric keyboard (T201). Once the wished value is set, press the **PRINT** button.

- It is now asked to weigh the glass float in air, confirm this with **PRINT**

dUE 1.81d

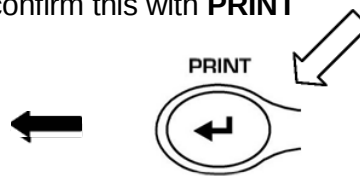


- Now tare if necessary and put the glass float to the weighing hook. Press then the **PRINT** button to acquire the value. During the acquisition of the value the message

perform a

WEI AIR will be flashing.

- It is now asked to weigh the glass float immersed into the liquid. Perform a tare if necessary and immerse the glass float into the liquid. Press the **PRINT** button. During the acquisition of the value the message **WEI LIQ** will be flashing.

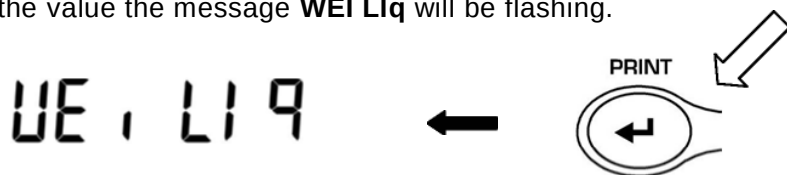

- The result of the density of the liquid is now displayed. Pressing the **PRINT** button it is possible to print the value of density, if the balance is equipped with a printer.



- In case of error the following message is displayed:




- Press the **ON/OFF** button to escape from density function, or the **MENU** button to perform another measurement.

20 Maximum load determination function

* The balance must be equipped with the specific optional software (non available for all models) *

The function **M LOAD** allows to measure the maximum load for a solid.

1. To activate the function **M LOAD** from zero display condition, press repeatedly the **MENU** button until the message “**M LoAd**” is displayed, then press the **PRINT** button to confirm. When this function is activated, a tare is automatically performed.

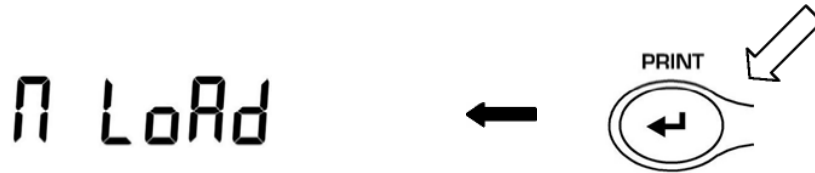

2. Activation of this function is indicated by the visualization of the letter **M** on the last digit on the display.



3. It is now possible to determine the weight of the maximum load.



4. Then press the **TARE** button to effect another measurement.



5. When the function **M LOAD** is active, the calibration is disabled.
6. Press the button **ON/OFF** to escape from **M LOAD** function.

21 RS232 interface features

1. General features

The balance transmits the value visualized on the display following serial RS232C standard, allowing to print the value of weight to a PC monitor or to a serial printer. In the case of connection to a PC, it will be possible to select the transmission in continuous mode or transmission at user command through pressing of the **PRINT** button (as described at pag.17). The balance is also capable of receiving commands, always through the standard RS232C, that allow to perform all the functions available through the keyboard of PC itself. The speed of transmission and reception can be selected, as described previously (pag.19), to 1200, 2400, 4800, e 9600 baud. The character format is of 8 bit preceded by one bit of start and followed by a bit of stop. Parity is not considered.

2. Selection of interface for PC

Selecting the transmission to PC (personal computer IBM compatible), it is achieved a continuous transmission output, at the same rate of weight update on display of the balance. It is possible to perform all the functions of the balances directly from the computer keyboard, transmitting to the balance the ASCII codes as in the table below. The connector to use for connection to PC is the number 1 (fig.1 pag. 37).


CODE	1 st FUNCTION (SINGLE PRESS)
"T" = H54	TARE
"C" = H43	CALIBRATION
"E" = H45	ENTER
"M" = H4D	MENU
"O" = H4F	ON/OFF

CODE	2 nd FUNCTION (PROLONGED PRESS)
"t" = H74	TARE
"c" = H63	CALIBRATION
"e" = H65	ENTER
"m" = H6D	MENU
"o" = H6F	ON/OFF

3. Selecting the transmission to PC at user command, it is achieved a transmission output only when the **PRINT** button is pressed, also in this case it is possible to perform all the functions of the balance directly from the keyboard of the computer, sending to the balance the ASCII codes in the table above. The connector to use for connection to PC is the number 1 (fig.1 pag. 37).

4. Connection of the balance to PC

To receive/transmit data, connect the connector 1 (fig.1 pag. 37) of the balance to the serial port of the PC as shown below:

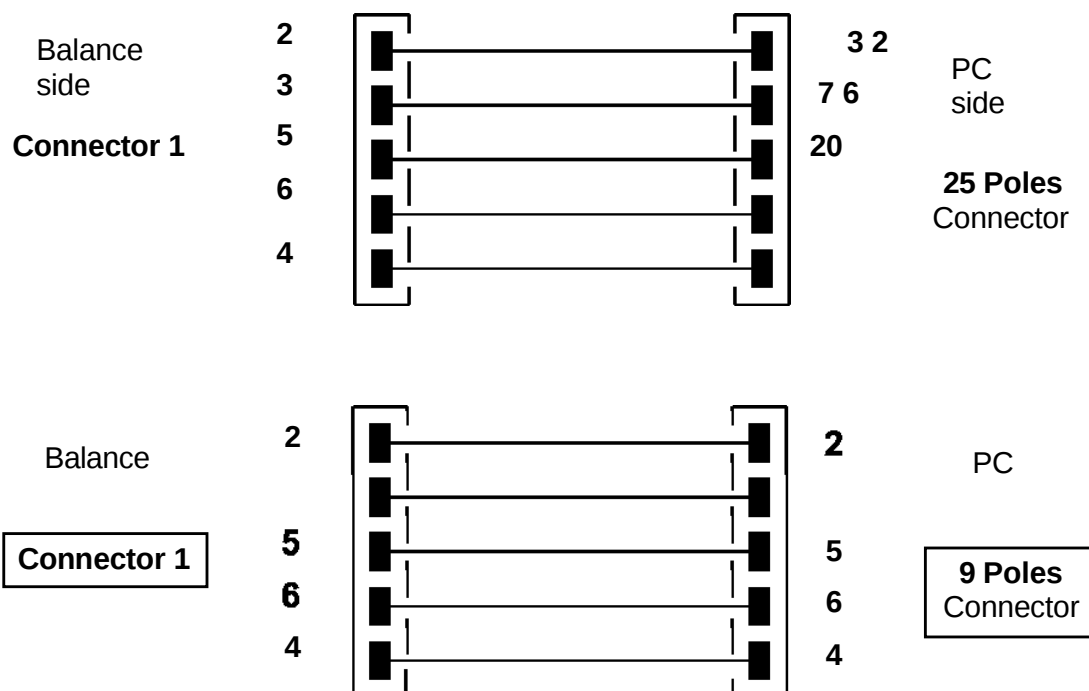

5. Transmission format

String transmitted is composed by the following 14 characters:

- ☐ First character: weight sign (blank or -)
- ☐ second/ninth character: weight or other data
- ☐ tenth/twelfth character: weight unit symbol
- ☐ thirteenth character: stability indicator
- ☐ fourteenth character: carriage return
- ☐ fifteenth character: line feed

Eventual non-significative zero are spaces.

6. In the following tables the various transmission formats are shown: Wei^ghin^g

mode (valid for continuous and at user command transmission modes)

1°	2°	3°	4°	5°	6°	7°	8°	9°	10°	11°	12°	13°	14°	15°r
Sign	Weight								Weight unit			Stability	CR	LF

Density mode (only in transmission at user command mode)

1°	2°	3°	4°	5°	6°	7°	8°	9°	10°	11°	12°	13°	14°	15°	16°	17°
d	=	Density value					Space		Weight unit						CR	LF

Piececounting mode (only in transmission at user command mode)

Number of pieces

1°	2°	3°	4°	5°	6°	7°	8°	9°	10°	11°	12°	13°	14°	15°	16°
Pcs			:	spaces				Number of pieces							

Total weigh of pieces

1°	2°	3°	4°	5°	6°	7°	8°	9°	10°	11°	12°	13°	14°	15°	16°	17°	18°	19°	20°
Weight						:	space	Value of weight								space	g	space	S

Unitary average weight of pieces

1°	2°	3°	4°	5°	6°	7°	8°	9°	10°	11°	12°	13°	14°	15°	16°	17°	18°
PMU			:	spaces				Value of weight							space	g	


7. Selection of interface to printer

Selecting the PRINTER mode, the serial output of the balance is set to work with serial printers.

In this case the printing is effected only when the **PRINT** button is pressed and with stable weight. If the stability is not reached within ten seconds, the message **ERROR05** is displayed preceded by a short acoustic signal and the value of the weight is not sent to the printer. The connector to be used for the connection is the number 1 (fig.1 pag.37).

8. Connection of the balance with the serial printer

Connect a serial printer to connector 1 (fig.1 pag.37) to the balance as shown in the following scheme:

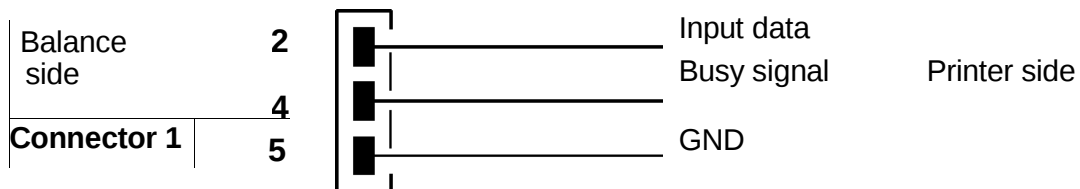

9. If the optional TLP50 printer is used, it will be possible to print both in *continuous-mode* and in *labels-mode* with the following formats:

Weighing mode and maximum load mode	Piececounting mode
--	---------------------------

12-02-2008 12:00 Weight: 22.000 g	12-02-2008 12:00 Pcs 100 Weight: 300.000 g PMU: 3.000 g
---	---

Density determination mode

12-02-2008 12:00 d= 2.80066 g/cm³d
--

10. Connection of the balance to the optional alphanumeric keyboard (code T201)

To connect the optional alphanumeric keyboard it must be used the connector 1 (fig.1 pag.37), the same used for connection to a PC. In this case the connection to PC or to printer must be effected through the connector placed on the optional alphanumeric keyboard.

22 Connectors places (rear)

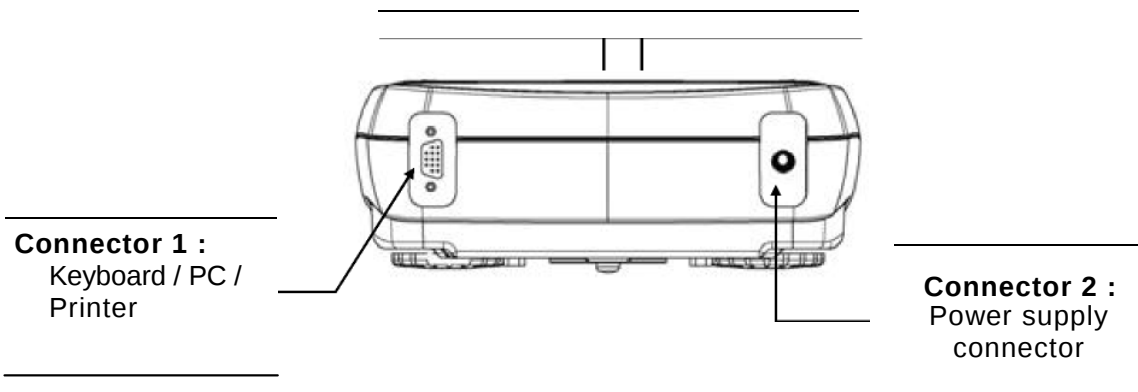


Fig. 1 Rear side of the balance

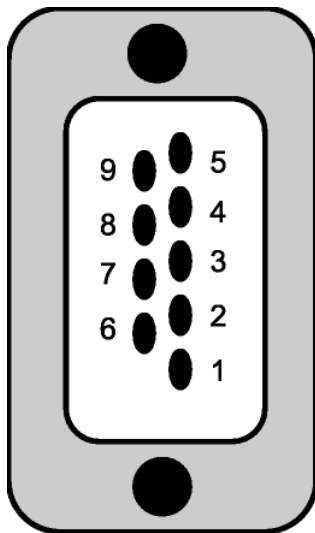


Fig. 2

CONNECTOR 1 PINS FOR KEYBOARD OR PC OR PRINTER

- ☐ pin 1 = Power +5V for keyboard
- ☐ pin 2 = Tx signal
- ☐ pin 3 = Rx signal
- ☐ pin 4 = busy signal
- ☐ pin 5 = Gnd
- ☐ pin 4-6 = connected to eachother for connection to PC

Fig. 3

23 Working with battery

Balance can work, at wish, using only the battery power.

The battery on board is located inside the balance and grants 10 hours of autonomy

1. With the external power supply disconnected, it is possible to switch on or switch off the balance trough the **ON/OFF** button and work without being connected to power AC outlet.
2. When the balance is working usingh the battery, the symbol of battery with charge level is displayed on up-right corner of the balance's display. The three lines all switched on means a maximum level of battery charge, instead when all three lines are switched off and only the battery symbol is displayed means that it is necessary to connect the balance to AC power outlet to re-charge the battery.



1. To re-charge the battery, connect the provided external power supply to the balance. (See Pag.5)
2. When the external power supply is connected, the balance automatically switch itself on and, after performing the initial test on display, it enters the stand-by mode showed by a single point in the display. In standby mode also the battery recharge symbol will remain switched on.



3. The charge level will step by step raise until it reaches the maximum level. However let the battery recharge at least for 8-10 hours before disconnecting the external power supply from the balance.


NOTE: Battery duration depends also from working mode of the display backlight. To get a longer battery duration, switch off the backlight or set it in AUTO mode.

24 Error codes

- ☐ **ERR01:** the weight does not reach stability after a tare operation ⇒ Protect the balance from air flows or from vibrations of the working table.
- ☐ **ERR02:** impossible to start the calibration operation due to balance instability ⇒ Protect the balance from air flows or from vibrations of the working table.
- ☐ **ERR03:** calibration weight not correct or balance unstable ⇒ Calibrate with correct weight or protect the balance from environment disturbs.
- ☐ **ERR04:** weight of samples for the piececounting function not adequate or unstable ⇒ Select a bigger number of samples or protect the balance from vibrations.
- ☐ **ERR05:** impossible to print due to instability ⇒ Protect the balance from environment disturbs.
- ☐ **ERR06:** the weight does not reach stability in density mode ⇒ Protect the balance from environment disturbs.
- ☐ **“UNLOAD”:** weight loaded on the pan or pan not positioned properly ⇒ Remove the weight from the pan or position properly the pan and underpan.
- ☐ **“CAL But”:** the balance requires to be re-calibrated ⇒ Unload weights, if any, on the pan, and press the CAL button.
- ☐ : Overrange condition ⇒ Unload the weights loaded on the pan.
- ☐ : Underrange condition ⇒ Place properly pan and underpan.


25 Quick guide to balance parameters setup

- To enter the balance parameters setup menu, press and keep pressed the **MENU** button until the acoustic alarm is over.
- Use then the **MENU** button to go to next parameterer, use the **CAL** button to go to previous and the **PRINT** button to cofirm the choice.
- To escape from menu, press the and keep pressed the **MENU** button until the acoustic alarm is over.

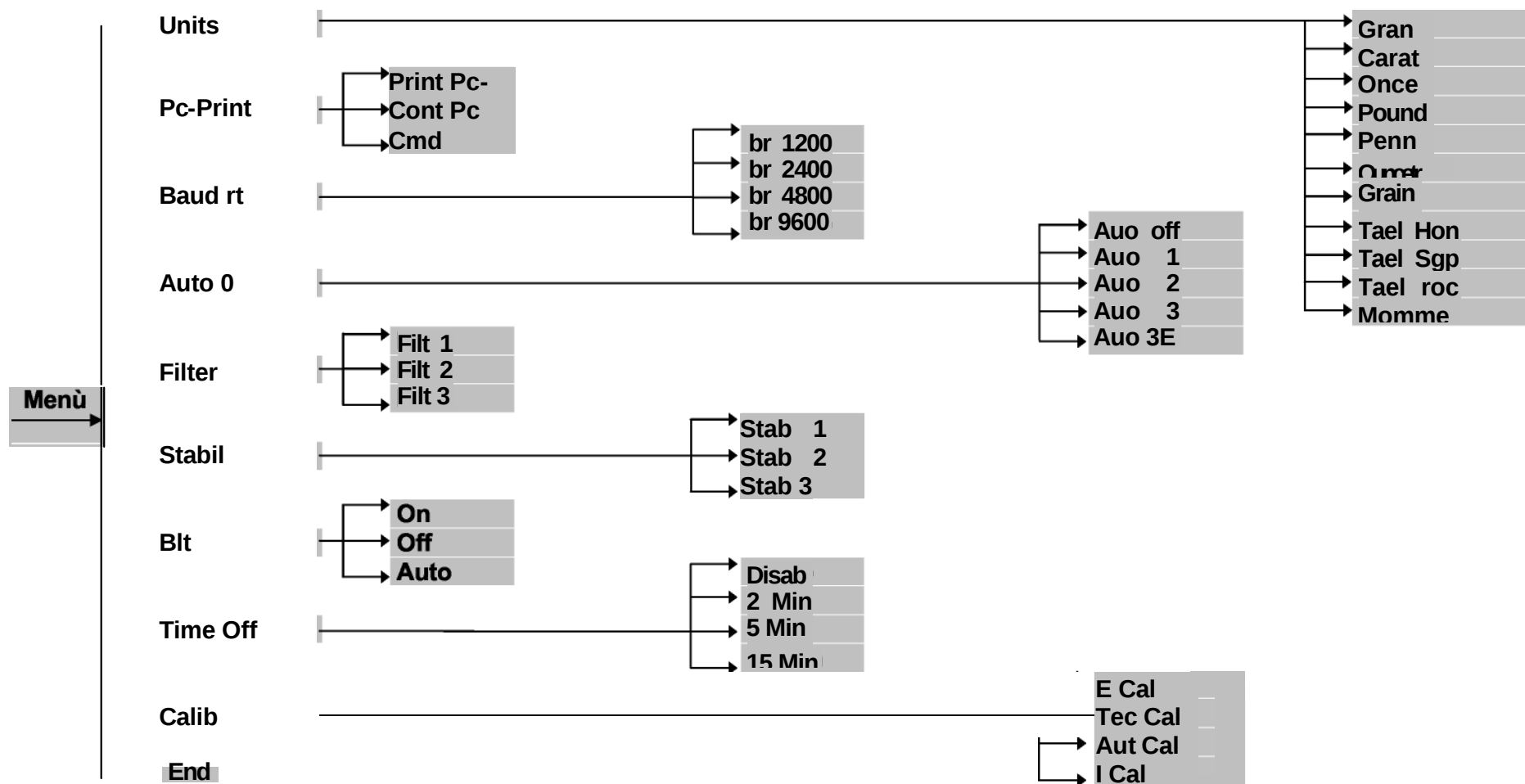

26 Quick guide to the use of the balance programs

- To enter the balance programs menu, press the **MENU** button.
- Use then the **MENU** button to go to next parameterer, use the **CAL** button to go to previous and the **PRINT** button to cofirm the choice.
- To escape from menu, press the and keep pressed the **MENU** button until the acoustic alarm is over.

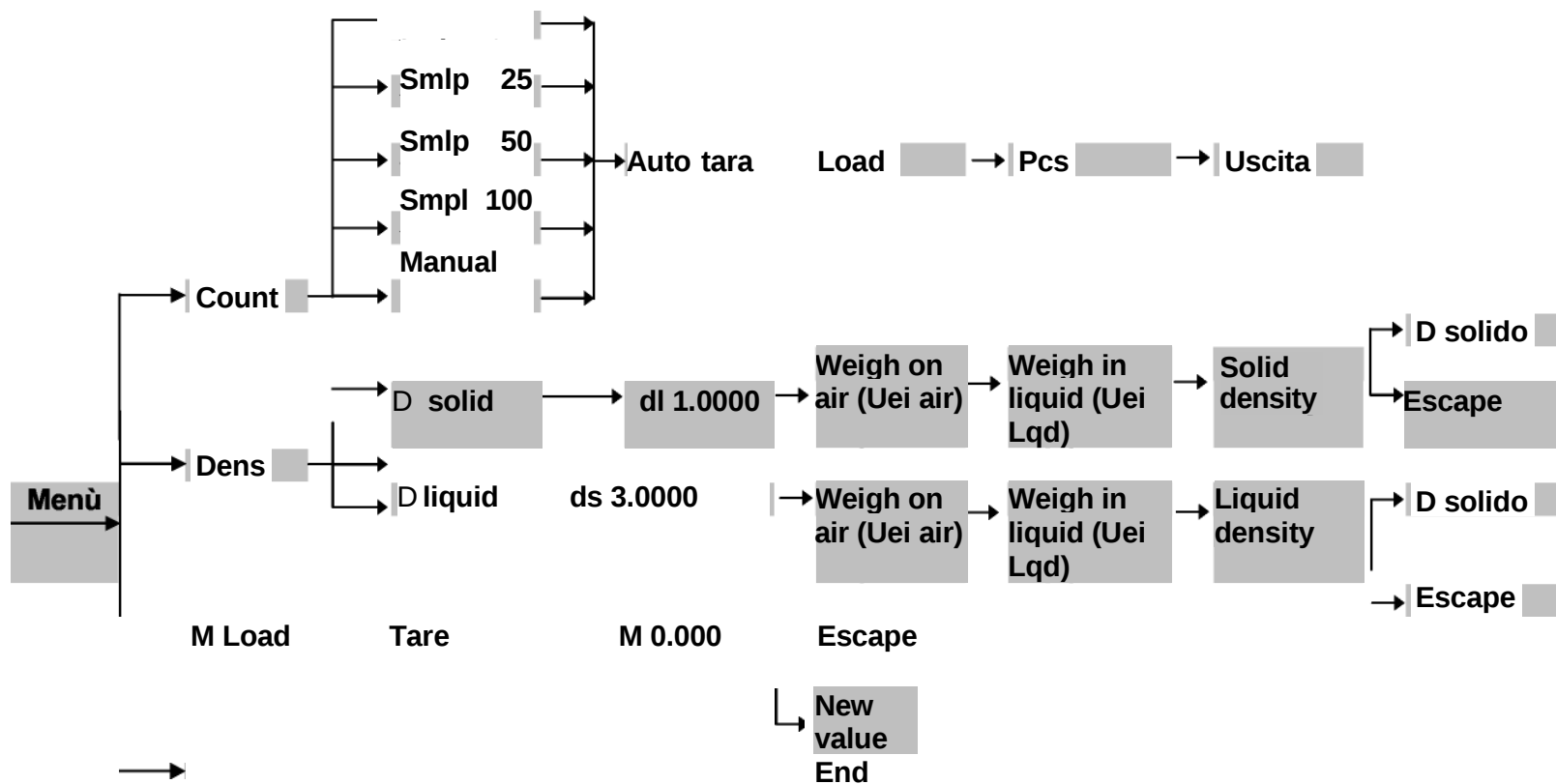

27 Warranty

- ☐ Duration of warranty is one (1) year from the date of purchase proved by invoice concerning the product or by delivery note.
- ☐ Warranty covers all parts resulting defective at the origin. It does not cover mechanical or electronic parts damaged by wrong installation, tampering or incorrect use.
- ☐ Warranty does not cover damages caused by impacts, balance drops or drop of objects on weighing pan.
- ☐ Shipment to and from service center is at customer charge

Intelligent Weighing Technology serves the industrial and laboratory markets in North America with competitive pricing and value for money on a complete range of industrial and laboratory scales and balances, force measurement and load monitoring technology from 0.01 mg to 275 tons.

- **We are the exclusive source for Intel I-Lab™, Intell-Industrial™, Intell-Check™, Intell-Count™, Intell-Weigh™, Intel I-Base™, Intel I-Scan™ and Intell-Print™ laboratory, industrial, counting, analytical, precision and toploading balances, printers, peripherals and weighing equipment.**
- **We are the exclusive source for UWE weighing equipment in the US and the Caribbean.**
- **We are the exclusive source for Intell-Lift™ Straightpoint (UK), Ltd. products in North America, Mexico and the Caribbean.**

We have a combined 50 years of experience in the weighing and measurement industry, both in the USA and worldwide. With contacts in over 50 countries, we provide you with the weighing and measurement equipment you need.

QUALITY- Products solidly built from the ground up with superior engineering and components for exacting results.

VALUE - Equipment priced for real-world business applications, with superior service and support.

EXPERIENCE - Expert advice to help you choose just the right product for your application.

Quality + Value + Experience...it adds up to the Intelligent Investment.

**Intelligent Weighing Technology,
Inc. www.intelligentwt.com**

