




GETTING STARTED WITH **TRACXER** **CODE READER**

TS201

PD801201

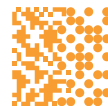

MICRONIC

Traceable Sampling Solutions

A CHECK COMPONENTS

Before installing the **Tracxer Code Reader TS201** be sure that you have all required items. Check the item list below!

- Tracxer Code Reader TS201;
- Single tube holder for 96-format tubes (already mounted);
- Serial-to-USB converter cable;
- CD-ROM with User Manual and Micronic Warranty Statement;
- This current document "Getting started".



If you do not have one of the items above, please contact the Micronic helpdesk by sending an email to: sales@micronic.com

B INSTALLATION

1 Installation conditions

Place the Tracxer Code Reader TS201 on a clean, dry, stable and horizontal surface.



Do not use the Tracxer Code Reader in an area with the following characteristics: high concentrations of dust, high humidity, explosion danger or toxic chemical storage.

2 Installation

- Connect the Tracxer Code Reader TS201 to the computer's USB port;
- Wait until the computer recognizes the reader;
- Open the program in which you want to use the scanned data.

C OPERATING TRACKXER CODE READER TS201

1 Operation

- Place a 2D Data-Matrix coded tube (any tube size in 96-, 48- or 24-well format) in the tube holder;
- The Tracxer Code Reader TS201 will scan and decode the tube code and export the code into the program which is activated on the computer (keyboard wedge functionality);
- The reader is set for robust reading of any 2D Data-Matrix code. An integrated light is switched on for easy and unambiguous sample identification when no external light is available.

2 Examples



Reads any tube size in 96-well format;



Reads any tube size in 48-well format;



Reads any tube size in 24-well format.

D WARRANTY

The warranty period on Micronic equipment begins on the day the product was purchased from an authorized dealer and ends after 12 months.

For more information about the Micronic Warranty on equipment please check the Micronic Warranty Statement file (PD801801) or paragraph 'Warranty and Liability' in the User Manual, which are both available on the supplied CD-ROM.



MICRONIC

Traceable Sampling Solutions



The working of the Traxcer Code Reader equipment is light sensitive. Despite of all precautions that have been taken by the designers of the system, there is always a possibility that exterior lighting conditions can interfere with the functioning of the Traxcer Code Reader. If the tube code is exposed too much to stray light, the correct functioning of the Traxcer Code Reader cannot be guaranteed.

PLEASE VISIT OUR WEBSITE: WWW.MICRONIC.COM

MICRONIC

Platinastraat 51
8211 AR Lelystad

Phone: +31 (0)320 277070
Fax: +31 (0)320 277055

E-mail: sales@micronic.com
Website: www.micronic.com