

WinScatter User Manual

Applied Photonics Group

Contents

1	Introduction	7
2	Installation procedures	9
3	Operating instructions	11
3.1	Connecting to the oscilloscope	11
3.2	Overview of the control panel and the main menu	12
3.3	Control panel	13
3.3.1	Scanning the waveform	13
3.3.2	Display options	14
3.3.3	Display operations	15
3.3.4	Cursors	15
3.3.5	Special functions	15
3.3.6	Data transfer	16
3.3.7	Waveform data	17
3.4	Main menu options	18
3.4.1	File	19
3.4.2	Edit	20
3.4.3	Operations	21
3.4.4	View	22
3.4.5	Help	22
3.5	Working off-line	22
4	Troubleshooting	25
4.1	Errors when running the application	25
4.2	Errors when using functions of the application	26
4.3	Errors in GPIB communications	27

List of Figures

3.1	Preparing the oscilloscope for external control through GPIB.	11
3.2	Connecting to the oscilloscope.	12
3.3	Control panel and main menu.	13
3.4	Data transfer between computer and oscilloscope.	16
3.5	Warning message.	17
3.6	Internal Memory Storage.	17
3.7	Special functions.	18
3.8	Parameters of the waveform.	18
3.9	Samples of the waveform.	19
3.10	Box displaying a waveform.	19
3.11	Main menu.	19
3.12	Dialog box for mathematical operations.	21
3.13	Explanation window	22
3.14	About window.	22
3.15	Off-line warning window.	23
3.16	Communication error message.	23

Chapter 1

Introduction

This manual is intended for operators who are responsible for operating the Sampling Optical Oscilloscope OOS-01 of Hamamatsu. This manual aims to:

- explain in detail how the application works,
- give the points to keep in mind when operating on the different devices for measuring the bandwidth in optical fibers, and
- advise how to obtain the highest performance.

This manual requires a user knowledge of Microsoft Windows operating systems as well as a working knowledge of terms related to optical fibers.

Furthermore, it is strongly recommended to read the user manual of the Optical Oscilloscope [1], since it is an extremely sensitive device and an incorrect usage could result in permanent damage.

Next, we are going to show the main features of the application and we will explain how it operates.

Chapter 2

Installation procedures

WinScatter works on Intel-based computers under Windows 95, Windows 98, Windows NT 4.0, Windows 2000 or Windows XP¹. The minimum requirements are the following: 16 MB RAM, 1.2 GB hard disk and a GPIB card.

In order to install the application, insert the WinScatter CD-ROM, select the WinScatter folder, and copy it into your computer. Next, open the WinScatter folder and run the batch file “RegisterMsFlxGrd.bat” in order to register the Microsoft FlexGrid ActiveX component (this action is absolutely necessary in order to be able to run the application). Afterwards, open the “WinScatter.ini” configuration file (using an ASCII text editor) and set the proper address parameters of your GPIB board and the Hamamatsu OOS-01 oscilloscope.

Once the application has been installed and properly configured, double click the “WinScatter.exe” application file and it will start up automatically.

¹The window style varies depending on the operating system. The figures shown later in this manual were taken using Windows XP.

Chapter 3

Operating instructions

3.1 Connecting to the oscilloscope

First of all, preparations have to be made in the oscilloscope so as to enable the external control through the GPIB interface. This is done following these steps:

1. Power on the oscilloscope and the laser driver.
2. Select CONT in the TRIG. MODE switch of the laser drive in order to input the trigger signal in the oscilloscope.¹
3. Using the control panel of the oscilloscope, press the SEL key twice and then the F4 (REMOTE) key as shown in Fig. 3.1.

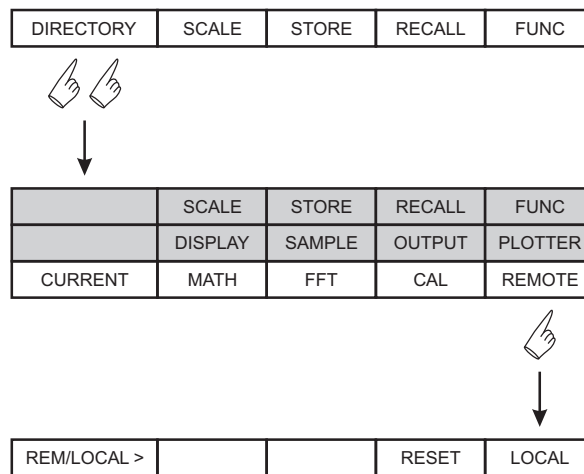


Figure 3.1: Preparing the oscilloscope for external control through GPIB.

Now the oscilloscope is in the remote control status.

If you now run the application, the window shown in Fig. 3.2 will be displayed. If the connection to the oscilloscope is successful, you will be able to choose between working on-line (all features enabled) or off-line (you can only access to certain functions). If the connection has failed, it can be due to one of the following three reasons:

¹From now on the laser diode is emitting. Therefore, special care should be taken.

- You are working on a computer which has no connection to the oscilloscope through GPIB. In that case, you will be able to work only off-line. It is important to note that this is not an error; this operation mode has been intentionally designed to allow the user to work with previously stored measurements in computers with no connection to the oscilloscope. This operation mode will be explained in detail in section 3.5 on page 22.
- The oscilloscope has not been correctly configured to enable external control through the GPIB card installed in the computer. Make sure that you have followed the instructions previously explained in this section. Then try to connect to the oscilloscope clicking the button “Try again”. After that, the application should display the button “On-line”.
- There is an error in the GPIB bus or the GPIB card installed in your computer. If you cannot connect to the oscilloscope in spite of having followed the steps explained before in this section, perhaps the GPIB cable is damaged or the GPIB card does not work properly. It is recommendable to note down the error code and refer to section 4.3 on page 27.

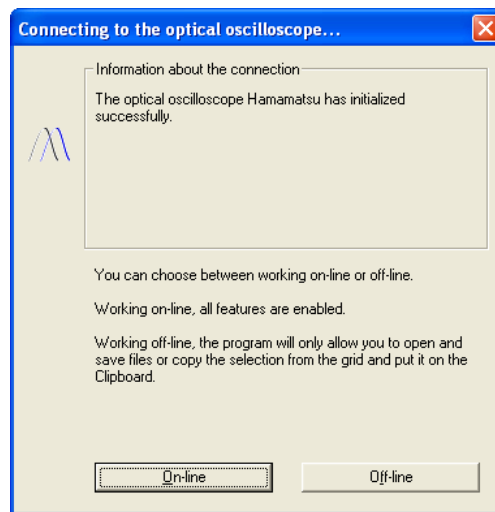


Figure 3.2: Connecting to the oscilloscope.

3.2 Overview of the control panel and the main menu

The control panel and the main menu are shown in Fig. 3.3.

The main options of the application are the following:

- **File.** It allows you to open or save files. You can also exit the application through this menu.
- **Edit.** It allows you to copy selected data from the grid “Waveform data” or export waveform data to other applications.
- **Operations.** This option allows you to display the dialog boxes corresponding to the data transfer between the computer and the oscilloscope, the internal memory storage

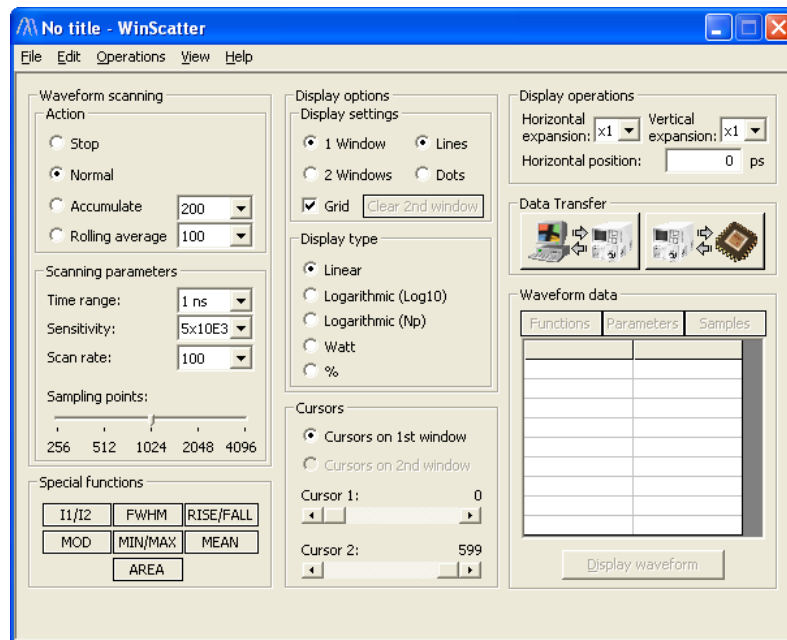


Figure 3.3: Control panel and main menu.

and the performance of mathematical operations with previously stored waveforms. It also allows you to switch off-line (the link to the oscilloscope is cut off).

- **View.** It allows to switch to the first or the second channel of the oscilloscope. It also allows to display the waveform obtained from the oscilloscope in the computer.
- **Help.** It displays a small window with operating instructions and gives information about version number and copyright.

Besides, the different options of the menu can be accessed quickly using the shortcut keys, such as, *Ctrl+Underlined Key*, *Alt+Underlined Key*, or the function keys (F1, F2, etc.).

3.3 Control panel

This section explains the operations available on the control panel and the menu.

3.3.1 Scanning the waveform

Inside the “Action” frame you can choose the scanning waveform method:

- **Stop.** It stops scanning waveform immediately.
- **Normal.** It starts scanning waveform.
- **Accumulate.** It accumulates waveform data. The combo box next to the radio button allows you to specify from 1 to 10000 accumulations.
- **Rolling average.** It averages waveforms while scanning. Each average is performed with the latest scan and the previously averaged results. The combo box next to the radio button allows you to specify from 1 to 100 averages.

The “Scanning parameters” frame contains the most important options that affect the way that waveforms are scanned. There are several combo boxes and sliders that allow customizing these parameters:

- **Time Range.** It sets the time range from 20 ps/DIV to 1 μ s/DIV.
- **Sensitivity.** Changes the sensitivity of the intensity scale. Depending on the type of display chosen (this issue will be explained below), the sensitivity is shown in watts or counts per division.
- **Scan range.** Sets the scan rate from 1 to 10000. If you increase the scan rate, there is an improvement in the signal-to-noise ratio, but special care must be taken, since: i) the acquisition time becomes longer (making the measurement results more sensitive to fluctuations in time and temperature), and ii) the internal circuitry of the photomultiplier tube is more sensitive to the light intensity (i.e. there is a decrease in the maximum value allowed for the light intensity incident on the oscilloscope). Therefore, if the message A/D SATURATED is displayed on the screen, the scan rate must be decreased.
- **Sampling points.** Sets the number of sampling points between 256 and 4096. The higher the number of sampling points, the higher the resolution of the waveform, but the higher the acquisition time.

3.3.2 Display options

The display settings allow you to modify how the waveforms are displayed on the screen of the oscilloscope. The options inside the frame “Display settings” are:

- **1 Window / 2 Windows.** It switches to the 1-window or 2-window displaying mode. In the former, the waveforms are displayed superimposed, whereas in the latter waveforms are displayed on separate windows.
- **Lines / Dots.** When dot-display is specified waveforms are displayed according to the samples acquired in the scanning process; whereas when line-display is specified, waveforms are displayed connected by lines.
- **Grid.** A grid showing the time and the intensity axes is displayed.
- **Clear 2nd window.** A click on this button clears the second waveform.

The radio buttons inside the frame “Display type” allow you to choose between a linear display and a logarithmic display:

- *Linear display:*
 - **Linear.** The waveform is displayed in a linear scale (the sensitivity is measured in counts per division).
 - **Watt.** The waveform is displayed in watts.
 - **%.** The waveform is normalized and it is displayed in %. The sensitivity cannot be changed.

- *Logarithmic display:*
 - **Logarithmic** (\log_{10}). The waveform is displayed in common logarithm (base-10 logarithm). The sensitivity cannot be changed.
 - **Logarithmic** (\ln). The waveform is displayed in natural logarithm. The sensitivity cannot be changed.

3.3.3 Display operations

The display operations allow you to modify the vertical and horizontal axes of the waveform, as well as to shift it along the horizontal time axis.

- **Horizontal expansion.** It expands the waveform in relation to the time axis. It can be expanded twice ($\times 2$), three times ($\times 3$) or four times ($\times 4$). The selection of the single expansion ($\times 1$) restores the original time axis.
- **Vertical expansion.** It expands or compresses the waveform relative to the intensity axis. This option cannot be modified if the logarithmic display mode is used.
- **Horizontal position.** You can shift the waveform display along the time axis by typing the time in picoseconds inside the edit box. With negative values the waveform shifts to the right, whereas with positive values it shifts to the left.

3.3.4 Cursors

The oscilloscope has two cursors that move along the time axis. The cursors are of great importance since they delimit the region where the special functions are applied. Each cursor is controlled by means of scroll bars placed at the bottom of the frame “Cursors”; the left side of the screen of the oscilloscope is specified by position 0, whereas the right side is specified by position 599.

Clicking the radio buttons placed above the scroll bars, you can choose between moving both cursors in the 1st window or in the 2nd window, provided that two waveforms are displayed.

3.3.5 Special functions

Clicking one of the buttons inside the frame “Special Functions” you can perform one of the seven special functions available. These functions are performed inside the limits defined by both cursors; therefore you should place correctly the cursors before performing these operations. The results are displayed on the grid located inside the frame “Waveform data” as explained in detail in subsection 3.3.7 on page 17.

- **I1/I2.** It determines the ratio between the intensities measured in the positions given by the respective cursors and the difference in time between them.
- **FWHM.** It calculates the Full Width at Half Maximum of a peak placed between both cursors.
- **RISE/FALL.** It determines the rise or fall time. The rise time is determined for the case when the intensity at the position of the cursor set at the left side of the screen is lower than the intensity at the position of the cursor set at the right side; otherwise, the fall time is determined.

- **MOD.** It determines the modulation factor of the area between cursors.
- **MIN/MAX.** It determines the maximum and minimum values in the area between cursors.
- **MEAN.** It determines the mean value of the intensities between cursors.
- **AREA.** It determines the accumulated value of the intensities between cursors.

3.3.6 Data transfer

The buttons inside the frame “Data transfer” allow you to exchange waveform data between the computer and the oscilloscope, or to store them in any of the four internal memories of the oscilloscope.

Data transfer between computer and oscilloscope

Figure 3.4 shows the button that pops up the “Data transfer” dialog box.

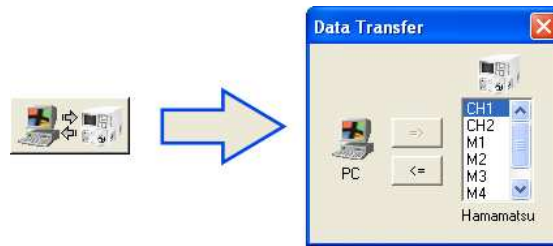


Figure 3.4: Data transfer between computer and oscilloscope.

By clicking the button that shows the arrow towards the right, the transfer takes place from the computer to the oscilloscope (upload), whereas if the button showing the arrow towards the left is clicked, the transfer takes place from the oscilloscope to the computer (download). There are six options available in the list box:

- **CH1.** When downloading, the waveform scanned by the photomultiplier tube connected to the port CH1 is transferred to the computer. When uploading, the waveform data is displayed on the 1st window.
- **CH2.** When downloading, the waveform scanned by the photomultiplier tube connected to the port CH2 is transferred to the computer. When uploading, the waveform data is displayed on the 2nd window.
- **M1, M2, M3, and M4.** The data transfer takes place between the computer and one of the four internal memories of the oscilloscope. If you request data from an empty memory, the application shows a warning message (see Fig. 3.5).

At the bottom of the control panel a progress bar shows the progress of the transfer. This dialog box can also be reached via the “Operations” menu.



Figure 3.5: Warning message.

Internal Memory Storage

Figure 3.6 shows the button that pops up the “Internal Memory Storage” dialog box.

When clicking the button that shows the arrow towards the right, the field “1st window” of the list box “Display Window” refers to the photomultiplier tube connected to the port CH1 of the oscilloscope, whereas the field “2nd window” refers to the tube connected to CH2. Therefore, the data storage takes place from one of the channels of the oscilloscope to the chosen internal memory.

Instead, when clicking on the button that shows the arrow towards the left, the data exchange takes place from the chosen internal memory to the chosen display window.

This dialog box can also be reached via the “Operations” menu.

3.3.7 Waveform data

The information corresponding to the waveform transferred to the computer is shown in the grid placed inside the frame “Waveform data”. The three buttons above the grid change the information type shown in the grid.

- **Functions.** Clicking this button the results for the special functions performed previously are shown. “No data” is displayed whenever a function has not been performed. An example is shown in Fig. 3.7.
- **Parameters.** Clicking this button the nine parameters of the waveform are shown (see Fig. 3.8 on page 18):
 - *Source.* It shows the source of the measured waveform sent to the computer: CH1, CH2, M1, M2, M3, or M4.
 - *Time Range.* It shows the time range used during waveform acquisition.
 - *Horizontal exp.* It shows the horizontal expansion used to acquire waveform data.
 - *Aperture time.* This parameter only makes sense when using CCD sensors instead of a photomultiplier tube.


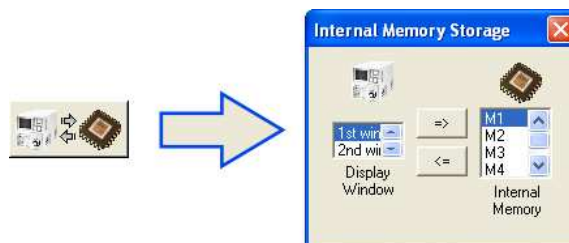


Figure 3.6: Internal Memory Storage.

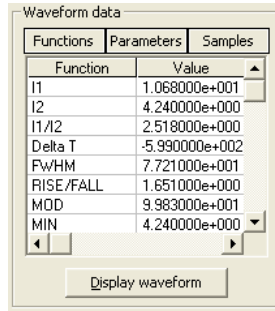


Functions	Parameters	Samples
Function	Value	
I1	1.068000e+001	
I2	4.240000e+000	
I1/I2	2.518000e+000	
Delta T	5.990000e+002	
FWHM	7.721000e+001	
RISE/FALL	1.651000e+000	
MDD	9.983000e+001	
MIN	4.240000e+000	

Display waveform

Figure 3.7: Special functions.

- *Sampling points*. It shows the resolution used during waveform acquisition.
- *Sensitivity*. It shows the sensitivity used to acquire waveform data.
- *Scan rate*. It shows the scan rate used during waveform acquisition.
- *Accum. count* It shows the total number of accumulated waveforms.
- *Channel no.* It shows the number of the photomultiplier tube used to scan the waveform. If there is only one tube connected to CH1, this parameter is always set to 0.

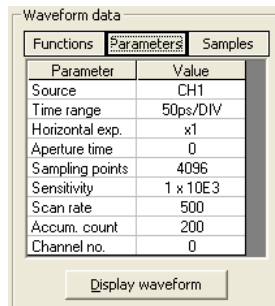

Functions	Parameters	Samples
Parameter	Value	
Source	CH1	
Time range	50ps/DIV	
Horizontal exp.	x1	
Aperture time	0	
Sampling points	4096	
Sensitivity	1 x 10E3	
Scan rate	500	
Accum. count	200	
Channel no.	0	

Display waveform


Figure 3.8: Parameters of the waveform.

- **Samples**. Clicking this button the samples of the measured waveform are shown. Each sample consists of a pair of values: the first value shows the time in ps and the second value shows raw data related to the number of intensity counts (these raw data do not take into account the selected scan rate and the number of accumulations). An example is shown in Fig. 3.9.

Finally, clicking the button “Display waveform” placed at the bottom of the grid, a box displaying the acquired waveform is shown (normalized to its maximum value). Figure 3.10 on page 19 shows the latter.

3.4 Main menu options

This section explains the options available through the main menu (those options already explained in the previous section will be briefly mentioned). Figure 3.11 shows the main menu.

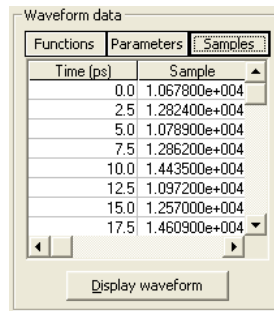

Time (ps)	Sample
0.0	1.067800e+004
2.5	1.282400e+004
5.0	1.078900e+004
7.5	1.286200e+004
10.0	1.443500e+004
12.5	1.097200e+004
15.0	1.257000e+004
17.5	1.460900e+004

Figure 3.9: Samples of the waveform.

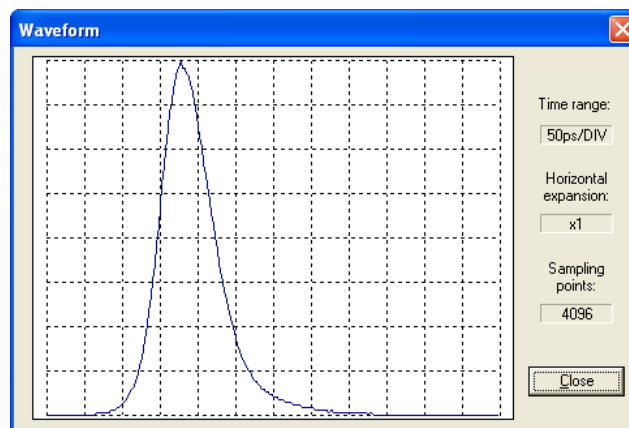


Figure 3.10: Box displaying a waveform.

3.4.1 File

It allows you to open or save files. You can also exit the application through this menu.

Open

It opens waveform data stored in a file with the ***.sct** default extension. If you attempt to open a file which does not contain valid data an error message will be displayed.

You can also access the dialog box “Open” using the shortcut key *Ctrl+O*.

Save and Save as...

The application stores the waveform data with the ***.sct** default extension. Selecting *Save*, you will overwrite previously saved data with the latest parameters, whereas selecting *Save*

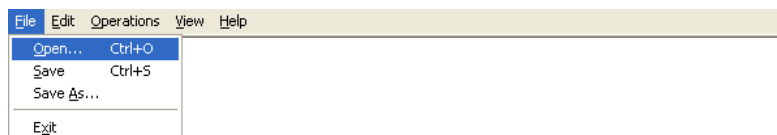


Figure 3.11: Main menu.

as... , the application will ask you a new name for the file to be created.

You can also access the dialog box “Save” using the shortcut key *Ctrl+S*.

Exit

You can exit the application in one of the following ways.

- Selecting “Exit” through “File” menu.
- Clicking the close special button on the right top of the window.
- Using the shortcut key *Alt+F4*.
- While maintaining *Alt* key, press *F*, release it and then press *X*.

To restore the control in the oscilloscope and, therefore, to make it possible to operate its front panel, you must press the F3 (RESET) key and then the F4 (REMOTE) key.

3.4.2 Edit

It allows you to copy selected data from the grid “Waveform data” or export waveform data to other applications.

It is also possible to display a pop-up menu with the same content as the “Edit” option with a right mouse click.

Export waveform to...

This option allows you to export the waveform data stored on your computer to other applications. WinScatter will automatically format the data in order to make them suitable for the application you have chosen to export to, and it will put them on the clipboard. In the following, there will be explained what you have to do to paste successfully the data on each application.

- **Spreadsheets.** WinScatter allows you to export waveform data to Microsoft Excel or OriginPro:
 - Microsoft Excel: open Microsoft Excel; create a new sheet, select the first cell and then click the button “Paste”.
 - OriginPro: open OriginPro; create a new workbook, select the first cell and then click the button “Paste”.
- **Mathematical applications.** WinScatter creates automatically the required code in order to export this to Mathematica (version 3.0 or higher) or MATLAB (version 5.1 or higher).
 - *MATLAB*. Open MATLAB; Select “M-file” through “File/New” menu and then the M-file editor will be displayed. Paste the content of the clipboard and change the name “winscatt” next to the *function* command with that one you will use to save the file. You can finally run it typing the file name in “MATLAB Command Window”.
 - *Mathematica*. Open Mathematica; paste the content of the clipboard on the command window and press “Intro”. Mathematica will automatically display the waveform.

Copy from grid

This option allows you to copy the selected data on the grid placed inside the frame “Waveform data” and put them on the clipboard. Then, they can be pasted in any other application, such as word processors or spreadsheets.

You can also copy the content using the shortcut key *Ctrl+C*.

3.4.3 Operations

It allows you to display the dialog boxes corresponding to the data transfer between the computer and the oscilloscope, the internal memory storage, and the performance of mathematical operations with previously stored waveforms. It also allows you to switch off-line (interrupting the communication with the oscilloscope).

Mathematical operations

If you select this option the dialog box shown in Fig. 3.12 is displayed.

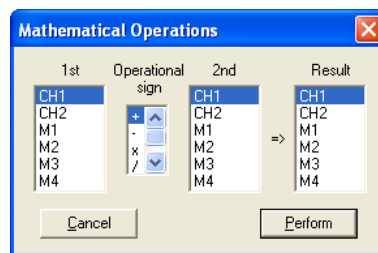


Figure 3.12: Dialog box for mathematical operations.

You can perform 4 different arithmetic operations on two waveforms. The measurement parameters (time range, sensitivity, number of samples, etc.) for both waveforms must be the same.

First of all, you have to select the first operand, then the operand (sum, rest, multiplication or division), then the second operand and finally the memory or window you want to store the result.

Clicking the button “Perform”, the oscilloscope performs the arithmetic operation. If you want to exit this dialog box without performing any operation, just click the button “Cancel”.

Data transfer

This is the same option as explained in subsection 3.3.6 on page 16.

Internal memory storage

This is the same option as explained in subsection 3.3.6 on page 17.

Off-line

Finally, this option allows you to switch off-line (the link to the oscilloscope is cut off). The latter will be discussed in section 3.5 on page 22.

3.4.4 View

This option allows to switch to the first or the second channel of the oscilloscope. It also allows to display the waveform obtained from the oscilloscope in the computer (the latter has been explained in subsection 3.3.7 on page 17).

3.4.5 Help

This option displays a small window with operating instructions and gives information about version number and copyright.

Figure 3.13 shows the operating instructions displayed when selecting “Operations”.

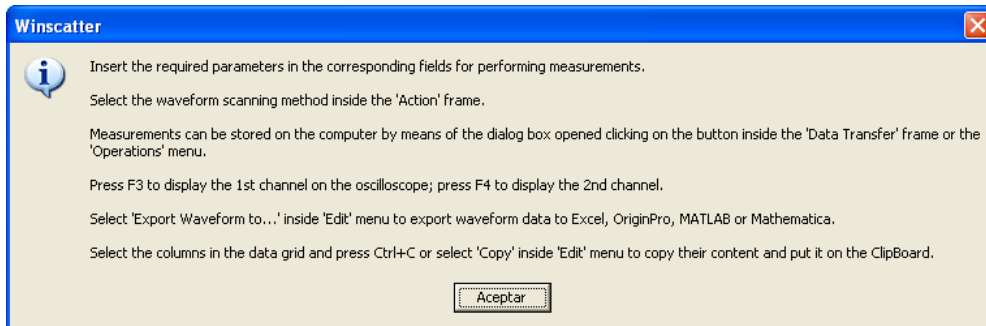


Figure 3.13: Window showing operating instructions.

Finally, Fig. 3.14 shows the window popped up when selecting “About WinScatter...”.

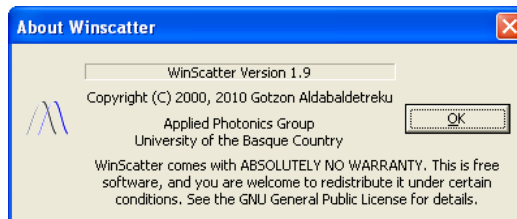


Figure 3.14: About window.

3.5 Working off-line

There are two ways to work off-line:

1. Clicking the button “Off-line” in the dialog box *Connecting to the optical oscilloscope* when you initialize the application (see Fig. 3.2 on page 12).
2. Selecting the option “Off-line” through the *Mathematical operations* menu (see subsection 3.4.3 on page 21). If you do this, the window shown in Fig. 3.15 will pop up.

When you work off-line, you can only access to a limited set of options:

- **Main Menu.** You can open and save files, display the waveform, and export waveform data to other applications.

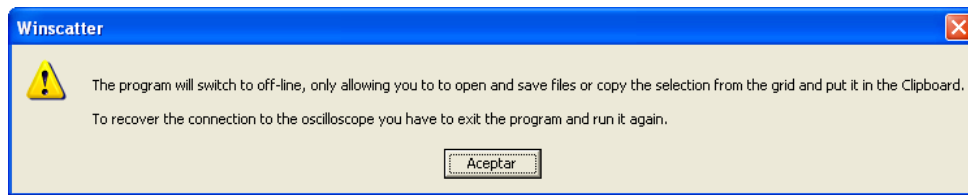


Figure 3.15: Off-line warning window.

- **Control panel.** You can only access to the controls located inside the frames “Special functions” and “Waveform data”. Working off-line, the special function buttons behave in the same way as the button “Functions” inside the frame “Waveform data”.

It is important to notice that when the communication link between the computer and the oscilloscope has been cut off, the application switches automatically to off-line, displaying an error message with the reasons for the failure (see Fig. 3.16).

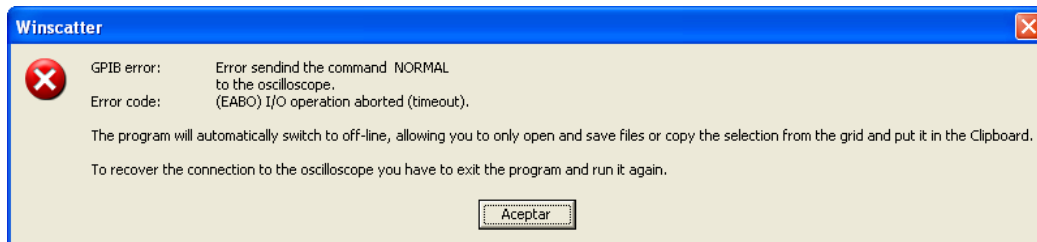


Figure 3.16: Communication error message.

In such a case, it is recommendable to note down the error code and refer to section 4.3 on page 27.

Chapter 4

Troubleshooting

4.1 Errors when running the application

- *The application does not run or there is a warning message showing there are not enough resources available.*

This means that there is not enough memory. The minimum required memory is 16MB.

If your computer fulfills these requirements but this message still appears, make sure there are no other applications running at the same time. Make also sure that there are at least 20MB free on the hard disk.

Although WinScatter works on the different 32-bit Windows operating systems, it is strongly recommended to use it on those systems based on NT technology (namely, Windows NT 4.0, Windows 2000 and Windows XP)¹.

- *When running the application there is a message in the “Connecting to the oscilloscope...” window showing that the connection has failed.*

This can be due to one of the following three reasons:

1. You are working on a computer which has no connection to the oscilloscope through GPIB. In that case, you can only work off-line. You should have a GPIB card installed on your computer and a GPIB cable connected to the oscilloscope.
2. The oscilloscope has not been correctly configured to enable external control through the GPIB card installed in the computer. Make sure that you have followed the instructions explained in section 3.1 on page 11. Then try to connect to the oscilloscope clicking the button “Try again”. After that, the application should display the button “On-line”.
3. There is an error in the GPIB bus or in the GPIB card installed in your computer. If you cannot connect to the oscilloscope in spite of having followed the steps explained before in this section, perhaps the GPIB cable is damaged or the GPIB card does not work properly. It is recommendable to note down the error code and refer to section 4.3 on page 27.

¹WinScatter has not been tested on Windows Vista or Windows 7 operating systems.

4.2 Errors when using functions of the application

- *When selecting any of the options of the control panel there is a message showing a GPIB error.*

This can be due to the following reasons:

1. You have gone back to the *LOCAL* mode pressing the *F4* key on the front panel of the oscilloscope, disabling the external control of the latter through the GPIB interface. For security reasons, the application will switch to off-line. If you want to restore the connection to the oscilloscope, you have to exit the application and run it again, preparing the oscilloscope by following the instructions explained in section 3.1 on page 11.
2. There is an error in GPIB bus or the GPIB card installed in your computer. Perhaps the GPIB cable is damaged or the GPIB card does not work properly. It is recommendable to note down the error code and refer to section 4.3 on page 27. In this case, the application will also switch off-line.

- *Sensitivity cannot be changed.*

The combo box “Sensitivity” is disabled when:

- The waveform scanning has been stopped with the command “Stop”.
- The execution of the commands “Accumulate” or “Rolling average” has finished.
- The “%” linear display type has been selected.
- Waveform data have been transferred from the computer to one of the channels (CH1 or CH2).
- The contents of one of the four internal memories has been transferred to one of the channels.

- *Vertical expansion cannot be changed.*

The combo box “Vertical expansion” is disabled when:

- There has been selected the “%” linear display type.
- There has been selected one of the two logarithmic display types *Logarithmic* (\log_{10}) or *Logarithmic* (\ln).

- *Horizontal expansion cannot be changed.*

The combo box “Horizontal expansion” is disabled when:

- The waveform scanning has been stopped with the command “Stop”.
- The execution of the commands “Accumulate” or “Rolling average” has finished.
- Waveform data have been transferred from the computer to one of the channels (CH1 or CH2).
- The contents of one of the four internal memories has been transferred to one of the channels.

- *The cursors cannot be changed to the 2nd window.*

In order to be able to select this option, it is necessary to have waveform data on the 2nd window. To achieve this you have to transfer data from the computer (see subsection 3.3.6 on page 16) or use the dialog box “Internal Memory Storage” (see subsection 3.3.6 on page 17).

- *The frame “Waveform data” does not display any information about the waveform displayed in the screen of the oscilloscope.*

It is necessary to transfer this waveform data from the oscilloscope to the computer (see subsection 3.3.6 on page 16).

- *When trying to transfer waveform data from one of the four internal memories to the computer there is a message showing “No data”.*

If you want to transfer data from one of the four internal memories it is necessary to have previously stored waveform data on that memory, transferring data from the computer (see subsection 3.3.6 on page 16) or using the dialog box “Internal Memory Storage” (see subsection 3.3.6 on page 17).

- *Cannot export waveform data.*

If you want to export waveform data to different applications, it is necessary to have waveform data on the computer. This can be achieved by transferring the waveform data from the oscilloscope to the computer (see subsection 3.3.6 on page 16) or by opening a file.

- *Cannot open a file or an error message is displayed.*

Since the application stores the waveform data with the `*.sct` default extension, you should only open this type of files (note that the application lets you open files with any other extension by selecting *All files (*.*)* in the list box). Opening a file which does not contain valid data will lead to displaying an error message.

It could also happen that the file is corrupted because the storage system has been damaged or any inconsistency in the File System of the operative system.

4.3 Errors in GPIB communications

In the following section the possible GPIB errors are enumerated and described. For more information, please refer to [2].

- *(EDVR) System error.*

This error occurs when you try to access a board or device that is not installed or configured properly.

- *(ECIC) Function requires GPIB board to be CIC.*

This error occurs when a command is sent to the GPIB board and the latter is not configured as the CIC or there are conflicts because of having multiple controllers.

- *(ENOL) No Listeners on the GPIB.*

This error usually occurs when a writing operation is attempted with no Listeners addressed. For a device writing, ENOL indicates that the GPIB address configured for that device in the software does not match the GPIB address of any device connected to the bus, that the GPIB cable is not connected to the device, or that the device is not powered on.

- *(EADR) GPIB board not addressed correctly.*

This error occurs when the GPIB board is CIC and is not properly addressing itself before the read and write functions. This error is usually associated with board-level functions.

- *(EARG) Invalid argument to function call.*

This error results when an invalid argument is passed to a function call.

- *(ESAC) GPIB board not System Controller as required.*

This error results when the functions requiring that the GPIB board be the System Controller are called and the latter has not this capability.

- *(EABO) I/O operation aborted (timeout).*

This error indicates that an I/O operation has been cancelled, usually due to a timeout condition.

- *(ENEB) Nonexistent GPIB board.*

This error occurs when no GPIB board exists at the I/O address specified in the configuration file. This occurs when the board is not physically plugged into the system, the I/O address specified during the configuration does not match the actual board setting, or there is a system conflict with the base I/O address.

- *(EDMA) DMA error.*

This error occurs if a system DMA error is encountered when the GPIB software attempts to transfer data over the GPIB using DMA.

- *(EOIP) Asynchronous I/O in progress.*

This error occurs when an asynchronous I/O operation has not finished before some other call is made. During asynchronous I/O, you can only use `ibstop`, `ibwait`, and `ibonl`, or perform other non-GPIB operations.

- *(ECAP) No capability for operation.*

This error results when your GPIB board cannot carry out an operation or when a particular capability has been disabled in the software and a call that requires the capability is made.

- *(EFSO) File system error.*

This error results when an `ibrdf` or `ibwrtf` call encounters a problem performing a file operation. Specifically, this error indicates that the function is unable to open, create, seek, write, or close the file being accessed.

- *(EBUS) GPIB bus error.*

This error results when a certain GPIB bus error occurs during device functions. All device functions send command bytes to perform addressing and other bus managements. Devices are expected to accept those command bytes within the time limit specified by the default configuration or the `ibtmo` function. EBUS results if a timeout occurred while sending these command bytes.

- *(ESTB) Serial poll status byte queue overflow.*

This error is reported only by the `ibrsp` function. ESTB indicates that one or more serial poll status bytes received from automatic serial polls have been discarded because of a lack of storage space.

- *(ESRQ) SRQ stuck in ON position.*

This error can only be returned by a device-level `ibwait` call with RQS set in the mask. ESRQ indicates that a wait for RQS is not possible because the GPIB SRQ line is stuck on.

- *(ETAB) Table problem.*

This error occurs only during the `FindLstn` and `FindRQS` functions. ETAB indicates that there was some problem with a table used by these functions.

Bibliography

- [1] Hamamatsu Photonics K. K., *Sampling optical oscilloscope OOS-01*. [Online]. Available: <http://www.hamamatsu.com/>
- [2] *GPIB user manual for Win32*, National Instruments, Austin, Texas, June 1998.