CLS-ARC 150 M

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Vielen Dank

CLS-ARC 150 M

This high quality product has been designed and developed to require a minimum of maintenance. Please follow the guidelines below and read this manual in its entirety to ensure safe operation of the unit.

Warning

- Please always disconnect mains plug before cleaning the device.
- Storing and using the Cold Light Source only in dry rooms.
- The provided power cord has a grounded plug. Please always use Power cable with correctly earthed socket.

Safety Information

- To avoid eye damage, don't look directly into the light source when turned on.
- To avoid burns, please let the Lamp cool down before replacing a spare part.

Symbols

The following symbols are used in this manual.

CAUTION! This symbol indicates a potential risk and alerts you to proceed with caution

Intended Use

The high performance cold light source CLS-ARC 150 M is particularly suited for image processing (machine vision). Further applications are be found in laboratories, automation, measurement technique, and in various industrial and technological lighting requirements.

The light source impresses by a modern and state-of-the-art design, which stands out by its modularity and its flexibility.

The CLS-ARC 150 M is a high quality light source, which provides the user a perfect light quality for an attractive price.

Package Contents

Description	Quantity
Cold Light Source CLS-ARC 150 M	1
Power Cord	1
Empty Filter Holder	1
User Manual	1

Getting Started

The Cold Light Source is a plug & play device very easy to use.


Please check before use whether ...

- ... the mains plug is connected.
- ... the main switch at the back side is switched on.
(blue LED at the front must glow)
- ... the lamp switch at the front side is switched on.
- ... the brightness is set to the desired value.
- ... the areas for the air intake (front) and out take (back) are free.
- ... the location has a good ventilation in order to avoid a thermal overload of the system.

Description of Buttons and Switches


Front View (fig.1)

- a) **Lamp switch**
To turn on and off the lamp.
- b) **Mechanical dimming with iris diaphragm**
For continuous dimming of the light without change the colour temperature.
- c) **Light Guide Access**
Connection of the light guide with a standard diameter of 15mm.
- d) **Air intake slots**
Air cooling must be ensured at any time.
- e) **Screw caps**
- f) **Cover of housing**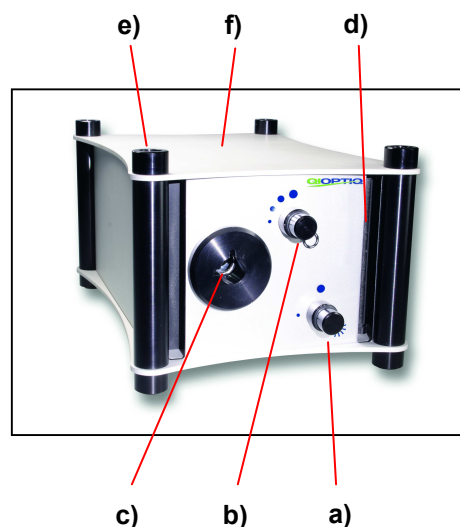


Back View (fig.2)

- g) **Plug for power cord**
- h) **Fine fuse**
The fine fuse can be removed by pulling out the insert with a small tool.
- i) **Main switch**
For switching on and off of all functions of the light source.
- j) **High quality low-noise rear fan**
The air out take must be ensured at any time.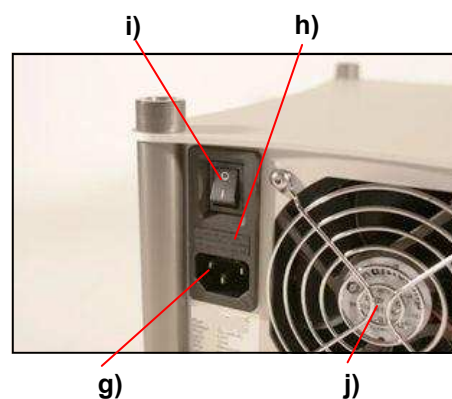

Interior (fig.3)

- k) **Cover plate**
- l) **Connecting clamp**
- m) **Lamp socket**
- n) **Halogen metal vapour lamp**
- o) **Lamp holder**

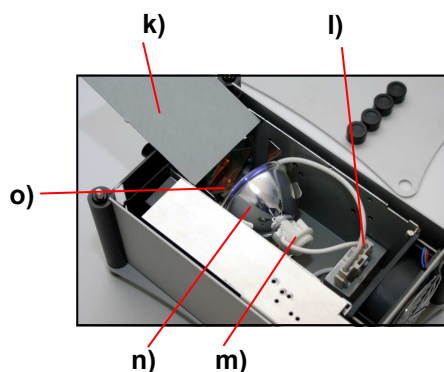CLS-ARC 150 M

Troubleshooting

The information in the table below can help to resolve the possible operating problems quickly.

Problem	Causes	Solution
No function at the device. No blue LED:	No network connection.	Please check, whether the power cord is connected.
	Fine fuse defective.	Please check the fine fuse (see picture). Replace the fuse in case of a defect. Please only use fuses with the right electrical values.
No light, but fan in operation. Blue LED is turned on:	Soiled air intakes could cause a too high temperature of the device. Automatic switch off by automatic thermo protection.	Please ensure a secure air intake. Clean the air intake and outtake areas.
	A too high environment temperature.	The environment temperature should not exceed 40° C max.
	Projection lamp is defective.	Please change lamp according to user manual.

Storage, Maintenance, Repair

The device is maintenance-free. Some points, however, you should take into consideration.

1. Please unplug the device before opening and cleaning.
2. For the cleaning of the housing please use a smooth, semi-wet cloth, if necessary with a smooth household cleaning fluid.
3. No moisture should entering into the interior of the housing.
4. Before replacing the spare lamp please clean the air intake and outtake areas with a small brush.
5. To avoid damages during the storing and transport please retain the packaging.
6. Repairs are only allowed by trained and experienced personnel. In case of in proper handling the warranty expires.

Technical Service

Web Resources:

Visit our Qioptiq website at www.qioptiq.com for:

- Contact details of technical customer service.
- Qioptiq product overview and price list as well as infos about accessories.
- Additional product information and special offers.

Contact:

For information or technical assistance contact directly Qioptiq Photinics GmbH & Co. KG.

Replacement of Lamps

1. Please switch off the halogen projection lamp with the electronic dimming switch (a).
2. Let the cold light source cool down for several minutes by using the fan.
3. Switch off the cold light source with the main switch (i).
4. Disconnect the power plug (g).
5. Unscrew the four screw caps at the surface of the device (e).
6. Remove the cover of the housing (f).
7. Open the cover plate (k).
Warning! The lamp could be hot. **Danger of burns!**
8. Ensure that the projection lamp has been cooled down.
9. Remove the projection lamp from the lamp holder (o).
10. Remove the lamp socket (m).
11. Please only use special halogen projection lamps 100W (Art.No.G40 5619 000) for replacement. **Attention!** The use of other lamps could cause damage in your device.
12. Clean the air intake and outtake areas (d and j) with a smooth tissue when replacing the lamp.
13. Please touch the new lamp only at the outside of the lamp.
14. Put the lamp socket (m) onto the new projection lamp.
15. Put the projection lamp in the lamp holder (o).
16. Close the inside cover plate (k).
17. Close the housing with the cover (f).
18. Fix the screw caps on the housing (e).
19. Plug in the power cord (g).
20. Now, your cold light source is ready to use.

Further repairs are only allowed by trained and experienced personnel. In case of improper handling the warranty expires.

Technical Specifications	
Weight:	3 kg
Dimensions:	260 x 175 x 135mm (L x W x H)
Operating voltage:	100-240 VAC / 50/60 Hz
Lamp:	150 W halogen metal vapour lamp
Average lamp life span:	6.000 h
Light guide access:	standard 15 mm
Light control:	Mechanic dimmer (iris diaphragm)
Color temperature:	4.200 K
Material:	Aluminium
Colour:	Silver, nature and titan anodized
Cooling system:	Active cooling by a low-noise-fan

CLS-ARC 150 M

Accessories											
	<table> <tr> <td>Daylight filter</td> <td>- Part. No.: G04 5610 000</td> </tr> <tr> <td>Green filter</td> <td>- Part. No.: G04 5611 000</td> </tr> <tr> <td>Yellow filter</td> <td>- Part. No.: G04 5612 000</td> </tr> <tr> <td>Red filter</td> <td>- Part. No.: G04 5613 000</td> </tr> <tr> <td>Blue filter</td> <td>- Part. No.: G04 5614 000</td> </tr> </table>	Daylight filter	- Part. No.: G04 5610 000	Green filter	- Part. No.: G04 5611 000	Yellow filter	- Part. No.: G04 5612 000	Red filter	- Part. No.: G04 5613 000	Blue filter	- Part. No.: G04 5614 000
Daylight filter	- Part. No.: G04 5610 000										
Green filter	- Part. No.: G04 5611 000										
Yellow filter	- Part. No.: G04 5612 000										
Red filter	- Part. No.: G04 5613 000										
Blue filter	- Part. No.: G04 5614 000										
	<p>Wide range of standard light guides and individual customize solutions.</p>										

Warranty

Qioptiq Photonics GmbH & Co. KG warrants that this product will be free from defects in material and workmanship for a period of one (1) year from date of purchase. If a defect is present, Qioptiq will, at its option, repair, replace, or refund the purchase price of this product at no charge to you, provided it is returned during the warranty period. This warranty does not apply if the product has been damaged by accident, abuse, misuse, or misapplication, or from ordinary wear and tear.

For your protection, items being returned must be insured against possible damage or loss. This warranty shall be limited to the replacement of defective products.

It is expressly agreed that this warranty will be in lieu of all warranties of fitness and in lieu of the warranty of merchantability.

Equipment Disposal

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste. Instead, it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

Thank you

Notizen/Notes