



# AS300 CAPTURE & PC ANALYSIS SOFTWARE MANUAL

ALARM  
MONITORING  
DATA LOGGING



AS300 Alarm and Monitoring Data Analysis - [04/03/2013 to 05/03/2013]

File Edit Reporting Window Help



Transmitter: 02

Serial Number: 2771

First Reading: 04/03/2013 20:21:10

Last Reading: 05/03/2013 20:00:45

No. of Readings: 95

Probe 1

Max Temp (°C): -74.4

Min Temp (°C): -77.7

Average Temp (°C): -76.4

Probe 2

Max Temp (°C): N/A

Min Temp (°C): N/A

Average Temp (°C): N/A

Filter


Start: 04 March 2013 20:21:10

End: 05 March 2013 20:00:45

Filter

Graph Scale break

Data Graph Alarms

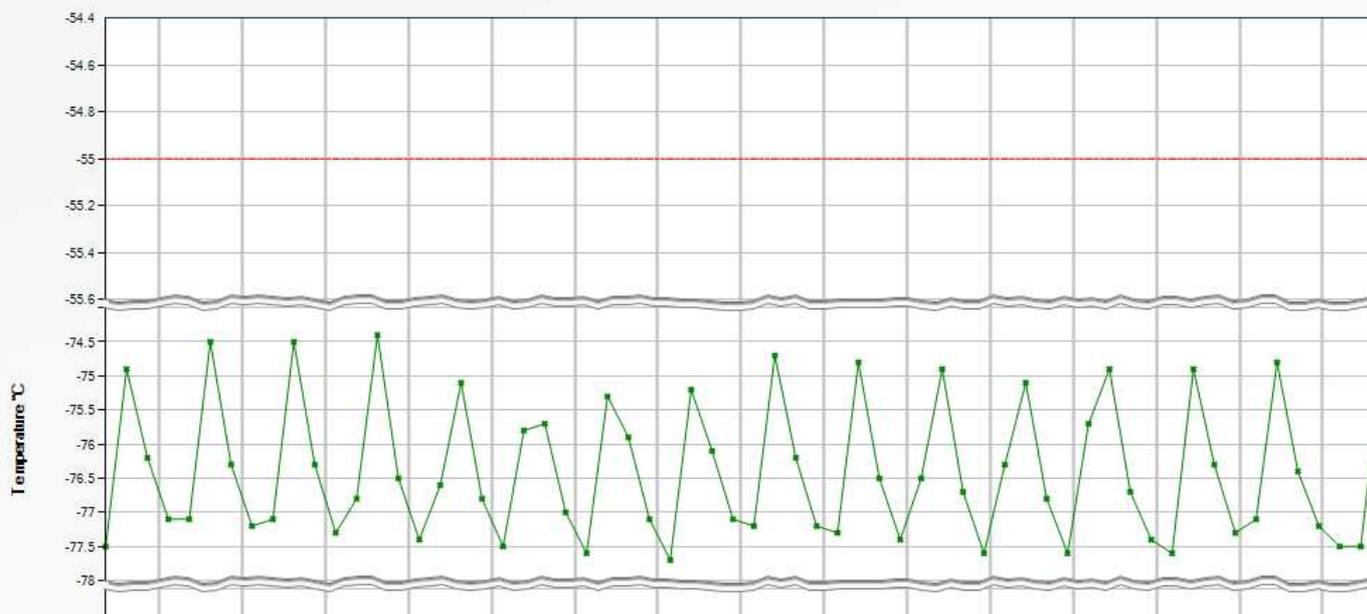

**Asper Systems Ltd**  
Laboratory Alarm & Monitoring Systems

## Table of Contents

**IMPORTANT:** THE LICENCE AGREEMENT SECTION 3 ON PAGE 7 MUST BE READ BEFORE INSTALLING OR USING SOFTWARE

1.0	Overview Capture Software	Page 3
1.1	Connecting a Computer to AS300	Page 3
1.2	Installing Capture Software	Page 3
1.3	Connecting to the AS300	Page 3
1.4	Main Screen	Page 3
1.5	Data Folders	Page 4
1.6	Descriptions	Page 4
1.7	Alarms	Page 4
2.0	Overview Data Analysis Software	Page 4
2.1	Installing Data Analysis Software	Page 4
2.2	Opening a Data File	Page 4
2.3	Using a Data File	Page 4
2.4	Top Screen Displayed Information	Page 4
2.5	Column Data	Page 4
2.6	Alarm Highlighted Column Data	Page 5
2.7	Filtering Data	Page 5
2.8	Displaying Graphs	Page 5
2.9	Graph Scale Break for Temperature	Page 5
2.10	X Axis Zooming	Page 5
2.11	Alarm Highlighting Graph Data	Page 6
2.12	Reporting	Page 6
2.13	Report Output Format	Page 6
2.14	Clipboard Export	Page 6
3.0	Licence Agreement	Page 7



## Section 1 ► Monitoring & Data Capture Software

### 1.0 Overview

- Display and logging of all the information sent from the transmitters.
- Data is updated every 15 minutes from each transmitter, increasing to every 5 minutes when in an alarm condition.
- Information sent from each transmitter (*If Enabled*) includes the following :- transmitter number, transmitter serial number, Tx Signal Strength, value I/P1, value I/P2, value I/P3, high alarm set-point, low alarm set-point, delay time and time of last update.
- Alarm status of input 1, input 2, input 3, AC power and low battery are displayed.
- Data is recorded in a CSV and PSA file. The PSA file is encrypted and can only be decoded and read by our analysis software. The CSV file can be read in our analysis software or excel. A set of new CSV and PSA files are generated every day to make record archiving simple.
- Standard 25 inputs can be expanded in blocks of 25 up to 100 inputs.
- Input I/P1, I/P2 and I/P3 descriptions can be added by customer.

Tx No	Serial No	Tx Signal	Description I/P 1	Description I/P 2	Description I/P 3	Value Input 1	Value Input 2	Value Input 3	High Alarm	Low Alarm	Delay Time	Alarm Condition	Last Update			
												STATUS	IP 1	IP 2	IP 3	AC
1												NO DATA				
2												NO DATA				
3												NO DATA				
4												NO DATA				
5												NO DATA				
6												NO DATA				
7												NO DATA				
8												NO DATA				
9												NO DATA				
10												NO DATA				
11												NO DATA				
12												NO DATA				
13												NO DATA				
14												NO DATA				
15												NO DATA				
16												NO DATA				
17												NO DATA				
18												NO DATA				
19												NO DATA				
20												NO DATA				
21												NO DATA				
22												NO DATA				
23												NO DATA				
24												NO DATA				
25												NO DATA				
26												NO DATA				

### 1.1 Connecting a Computer to AS300RX

Connect the supplied AS300 cable from the computer USB to the AS300. If drivers are required on your PC they can be downloaded from our web site at <http://www.asper.co.uk/Downloads.htm> **AS300 DRIVERS**, they are also supplied on the AS300 software pen drive stick with your Capture software.

### 1.2 Installing Capture Software

Double click on the Install file, prompts will guide you through the installation. Once installed the program can be accessed from the program bar under the heading Asper or screen icon.

### 1.3 Connecting to the AS300RX

The program is required to be run in the "administrator condition", to do this right click the mouse and click *Run as administrator*. Once the program has opened click on Settings > Comm port. Click the box Automatically Find USB Dongle, if the comm port is known this can be selected in Manually Select Comm Port. If more than 20 transmitter are on the system please Select Disconnected Detect Time to 30 mins or greater. Once this has been done go to the Settings > Connect to connect the AS300 to the PC. When connected the relevant Comm port will be displayed at the bottom RHS of the screen, on the LHS of the screen Asper Device Connected will be displayed. This procedure only needs to be done on the first start up of the Capture Software or if a different USB port is selected. With future restarts the capture software will connect automatically to the AS300

The Capture Software must be continually run to capture and store data .

If the program is not run in the Administrator Mode it may have problems with disconnections from the PC to the AS300.

### 1.4 Main Screen

The data will collect the following details from each transmitter, transmitter number, transmitter serial number, transmitter signal, value I/P1, value I/P2, value I/P3, high alarm set-point, low alarm set-point, delay time and time of last update. This data will be collected every 15 minutes, this interval will increase to every 5 minutes when the transmitter is in an alarm condition. If a probe is broken at a transmitter PB will be displayed at the input value.

### 1.5 Data Folders

The user can enter the path of two data folders where the collected data will be stored. File > Data Folders > Data Folder1 + Data Folder 2.

### 1.6 Descriptions

The user can enter descriptions for I/P1, I/P2 and I/P3 for each transmitter. Most transmitters purchased only use I/P1 and a typical description would be type of equipment or ID or location. E.g. -80 Freezer No 05.

### 1.7 Alarms

Alarms are indicated on the screen by the colour of the transmitter monitoring box, changing from green to red. Items monitored are I/P1, I/P2, I/P3, AC connected and battery low.

## Section 2 ► DATA ANALYSIS SOFTWARE

### 2.0 Data Analysis Software Overview

- CSV and encoded PSA files can be read.
- Numerical & Graph information easily displayed for any transmitter.
- Easy print facility for data required.
- Filtering of data is possible between specific dates and times.
- Statistical Information of the following is recorded : First Reading; Date and Time, Last Reading; Date and Time, Number of Readings, Maximum Temperature, Minimum Temperature, Average Temperature, Time in High Alarm and Time in Low Alarm.
- The Data Capture software and Data Analysis software come as a package. Each licence allows one capture software and up to 5 data analysis software programs.

### 2.1 Installing Data Analysis Software

Double click on the AS300 Alarm and Monitoring Data Analysis file, prompts will guide the installation. Once installed the program can be accessed from the program bar under the heading Asper.

### 2.2 Opening a Data File

To open a data file go to File > Open. The files are automatically saved in a date form e.g. 12\_05\_2013 is data collected on the 12th May 2013. Two files are made for each date a csv and psa. The psa file is encrypted and can only be viewed in the analysis software.

### 2.3 Using a Data File

Once the required date or dates are opened for analysis, select the transmitter unit number required to be viewed. This is done by using the mouse to move the "Transmitter Unit" column numbers up or down. Once the required transmitter unit is selected data for this unit is displayed.

### 2.4 Top Screen Displayed Information

Information at the top of the screen is displayed for the selected transmitter. The information is Serial Number, Descriptions if entered into the capture screen, First Reading, Last Reading, Number of Readings, Max Temp, Min Temp and Average Temp for I/P1 (Probe 1) and I/P2 (Probe 2)

### 2.5 Column Data

The data in the columns gives easy access to the following :-

Date, Time, Temperature input 1, Temperature Input 2, Door Switch I/P3, AC Power, Battery Condition, High Alarm Set-point, Low Alarm Set-point and Delay Time.



### 2.6 Alarm Highlighting Column Data

If an alarm occurs it will be highlighted in red for a high alarm and blue for a low alarm in the column data.

Date	Time	Temperature	High TCO	Low TCO	High Alarm	Low Alarm	High Alarm	Low Alarm
24/02/2007	08:04:41:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:05:01:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:24:48:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:24:49:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:24:50:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:24:51:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:24:52:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:24:53:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:24:54:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:24:55:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:24:56:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:24:57:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:24:58:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:24:59:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:00:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:01:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:02:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:03:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:04:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:05:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:06:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:07:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:08:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:09:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:10:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:11:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:12:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:13:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:14:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:15:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:16:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:17:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:18:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:19:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:20:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:21:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:22:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:23:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:24:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:25:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:26:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:27:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:28:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:29:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:30:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:31:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:32:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:33:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:34:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:35:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:36:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:37:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:38:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:39:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:40:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:41:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:42:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:43:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:44:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:45:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:46:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:47:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:48:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:49:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:50:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:51:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:52:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:53:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:54:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:55:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:56:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:57:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:58:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:30:59:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0
24/02/2007	08:31:00:012	0.0	0.0	0.0	0.0	0.0	0.0	0.0

### 2.7 Filtering Data


Data can be filtered by ticking the filter box and selecting a start date and time and end date and time. Once this has been done click on the filter button.

### 2.8 Displaying Graphs


A graph of the current data can be displayed by clicking on the Graph Tab. To return to the column data, click on the Data Tab.

### 2.9 Graph Scale Break for Temperature

When the graph is displayed, if there is a relatively large difference between the High Alarm, Low Alarm and I/P1, I/P2 measurements, the I/P1 and I/P2 variations can be hard to distinguish. To make this clearer, the "Axis Scale Break" feature collapses the gap between the highest data line and lowest data line. This allows the variations between the data lines to become more visible.

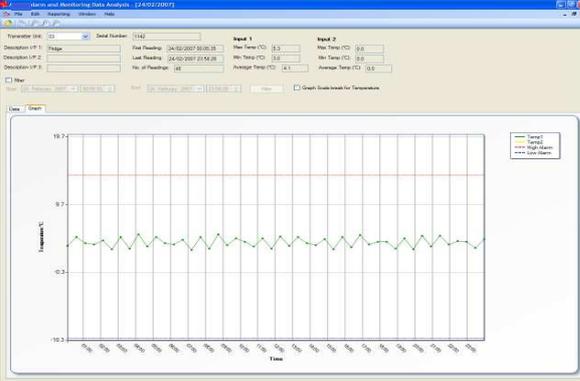


Scale Break Off

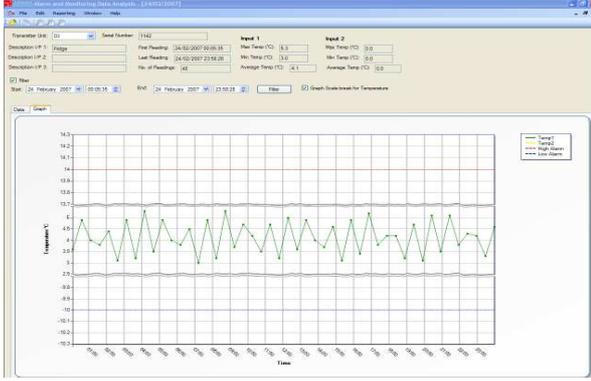


Scale Break On

### 2.10 X Axis Zooming

As well as the existing "Filter" functionality of the analysis application, which allows the user to filter a subset of the transmitter readings using a smaller time span, there is also the "Graph Zoom" function. This allows the user to zoom into a specific X-axis range of the graph. This is achieved by clicking on the start of the range required and then drag-

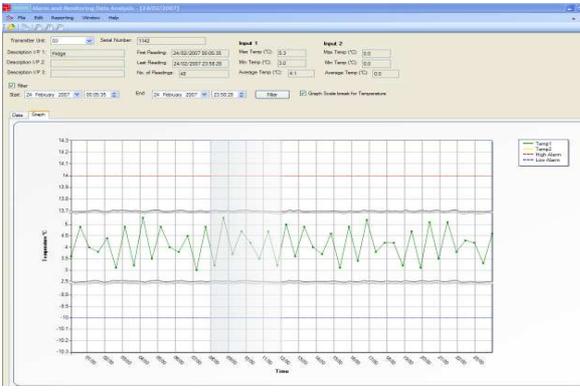


Highlighted Area to be Zoomed

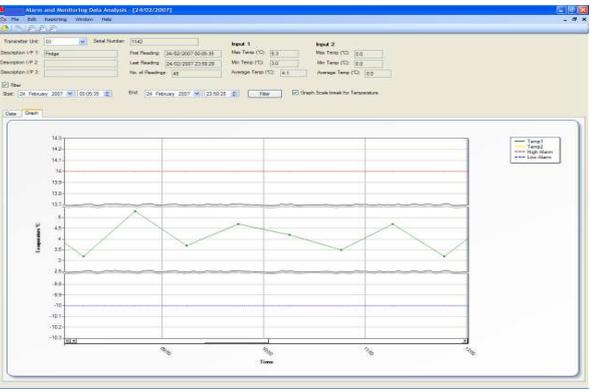

Zoomed Graph

ging the mouse to the end of the range the zoomed data is then displayed.

### 2.11 Alarm Highlighting

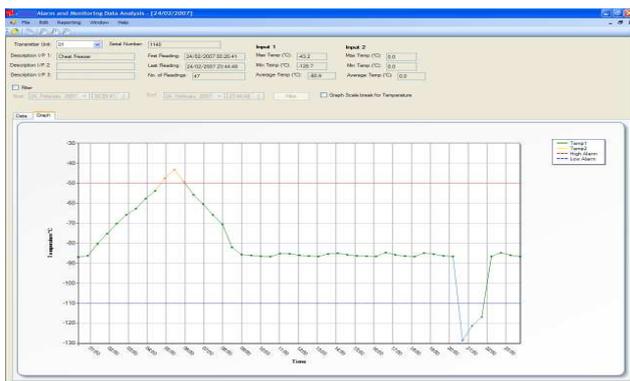

If either I/P1 (Temp1) or I/P2 (Temp2) value goes into the high alarm, the graph line is coloured red. If the input value goes below the low alarm set-point, the graph line is coloured blue.


### 2.12 Reporting

Print out of data can be done in three ways.

- “Transmitter Report” prints the data column view.
- “Graph Report” prints the graph view.
- “Summary Report” prints the “header” information for each transmitter.

To access these functions go to Reporting > Transmitter Report etc.

Transmitter Report

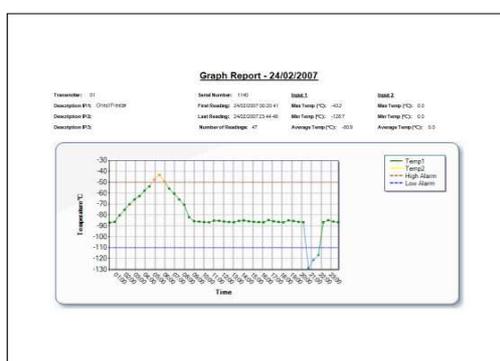

Graph Report

Summary Report

### 2.13 Report Output Format

The reports can be output in the following formats: -

- Adobe Acrobat (PDF)
- Microsoft Excel
- Microsoft Word



This can be done by selecting the “Exporting Report” button from the top left of the report window. Then select the report type from the “Save as type” selection in the “Export Report” dialog window.

### 2.14 Clipboard Export

Data in the data column and graph can be copied to the windows clipboard. In the data column the user can select the appropriate data cells using the mouse. The user then select the “Edit / Copy” menu item from the main window. For graphs the Graph’s image is copied to the Windows Clipboard  
 Once the data is copied to the Windows Clipboard, this can then be pasted into any appropriate Windows application  
 (Note :- By clicking the top left cell, all the cells in the data column are selected)

## Section 3 ► CAPTURE AND ANALYSIS SOFTWARE COPYRIGHT

Asper Capture and Analysis Software Copyright © 2013 by Asper Systems Ltd  
All rights Reserved

This license applies to the standard-licenced version of Asper Capture and Analysis Software.  
**Asper System Ltd Software Licence**

### Your Agreement to this Licence

You should carefully read the following terms and conditions before using, installing, copying, or distributing this software. Unless you have a different license agreement signed by Asper Systems Ltd, your use, installation, copying, or distribution of the Capture and Analysis Software indicates your acceptance of this agreement ("Licence"). If you do not agree to all of the terms and conditions of this Licence, then: (a) do not use, install, copy, or distribute any copy of the "Licence" Software with which this Licence is included.

### Scope of Licence

The terms and conditions of this Licence describe the permitted use and user's) of each Licenced Copy of Asper Capture and Analysis Software. Each licence allows the organization or user to one installation of the Capture Software and up to five installations of the Analysis Software.

All rights of any kind in the "Licenced" software which are not expressly granted in this Licence are entirely and exclusively reserved to and by Asper Systems Ltd. This software is protected by copyright laws and other applicable laws and treaties. You may use, install and distribute the "Licenced" software solely as expressly provided in this Licence. You may not rent, lease, loan, sublicense, modify, translate, reverse engineer, decompile, disassemble, or create derivative works based on, the Capture and Analysis software, nor permit anyone else to do so.

### Warranty Disclaimers and Liability Limitations

**Asper Capture and Analysis Software, and any and all accompanying software, files, data and materials, are distributed and provided "AS IS" and with no warranties of any kind, whether express or implied, including, without limitation, any warranty of merchantability or fitness for a particular purpose. Asper does not warrant, guarantee, or make any representations regarding the use of, or the results of the use of, The Capture and Analyst Software. Asper Systems Ltd does not warrant that the operation of the Capture or Analysis software will be uninterrupted or error-free, or that the use of any passwords and /or encryption features will be effective in preventing the unintentional disclosure of information contained in any file.** Further, there is no warranty for the optional virus scanning feature. (The Capture and analysis software does not scan for viruses; it simply runs external programs that claim to perform this function.) You acknowledge that good data processing procedure dictates that any program, including the Capture and Analysis Software, must be thoroughly tested with non-critical data before there is any reliance on it. **This disclaimer of warranty constitutes an essential part of this Licence.**

**Except to the extent expressly prohibited by applicable statutes, Asper Systems Ltd shall not be liable for any damage or financial loss arising out of the use of, or inability to use, the Capture and Analysis Software. Any liability of Asper Systems Ltd will be limited exclusively to refund of the licence fee paid to Asper Systems Ltd. Except to the extent expressly prohibited by applicable statutes, in no event shall Asper Systems Ltd, or its principals, shareholders, officers, employees, affiliates, contractors, subsidiaries, or parent organisations, be liable for any indirect, incidental, consequential, special, or punitive damages whatsoever relating to the use of the Capture or Analysis software, or to your relationship with Asper Systems Ltd, (including, without limitation, loss or disclosure of data or information, loss of profit, revenue, business opportunity or business advantage, or business interruption), whether based upon a claim or action of contract, warranty, negligence, strict liability, contribution, indemnity, or any other legal theory or cause of action, even if advised of the possibility of such damages. In addition, in no event does Asper Systems Ltd authorise you or anyone else to use the Capture and Analysis Software in applications or systems where the failure to perform can reasonably be expected to result in a significant physical injury, or in loss of property, or in loss of life. Any such use is entirely at your own risk, and you agree to hold Asper Systems Ltd harmless from any and all claims or losses relating to such unauthorized use.**

### General

This Licence is the complete statement of the agreement between the parties on the subject matter, and merges and supersedes all other or prior understandings, purchase orders, agreements, and arrangements. This Licence shall be governed by the laws of Scotland. Exclusive jurisdiction and venue for all matters relating to this Licence shall be in courts and for a located in Scotland, and you consent to such jurisdiction and venue. There are no third-party beneficiaries of any promises, obligations, or representations made by Asper Systems Ltd herein. Any waiver by Asper Systems Ltd of any violation of this Licence by you shall not constitute, nor contribute to, a waiver by Asper Systems Ltd of any other or future violation by you of the same provision, or any other provision, of this Licence. If any part of this Licence or the application thereof to any person or circumstance is for any reason held invalid or unenforceable, it shall be deemed severable, and the validity of the remainder of this Licence, or the applications of such provision to