


Operation Manual

Model VTX-3000L

Thank you very much for choosing the LMS Model VTX-3000L Mixer Uzusio.


Your Mixer Uzusio is CE Certified and designed with function, reliability, and safety in mind. For safe operation, please pay attention to the following:

- Check that the electrical supply from your power socket corresponds to the voltage stated on the Mixer Uzusio.
- Place your Mixer Uzusio on a clear, flat and level surface
- Never immerse your Mixer Uzusio in water; serious injury from shock or fire may result.

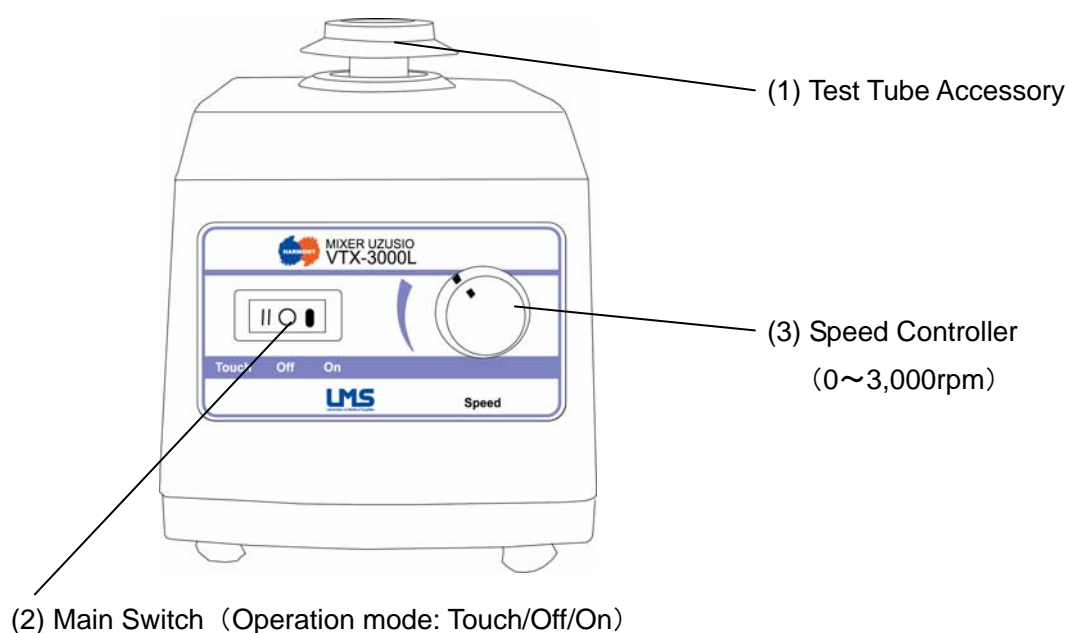


The unit ships with the following accessories.

- a. VTX-3000L Mixer Uzusio
- b. VTX-3000TT Test Tube accessory
- c. VTX-3000FP Flask accessory

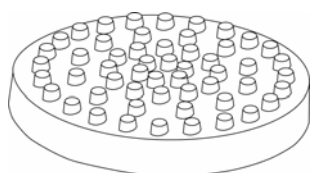


Optional Plates

- a. VTX-3000MP Microtitre plate Accessory

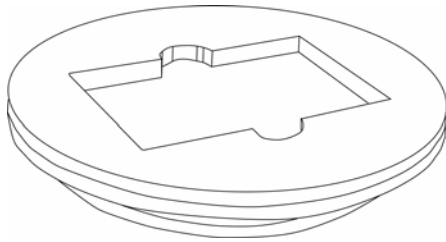

- b. VTX-3000MT Micro Centrifuge tube Accessory

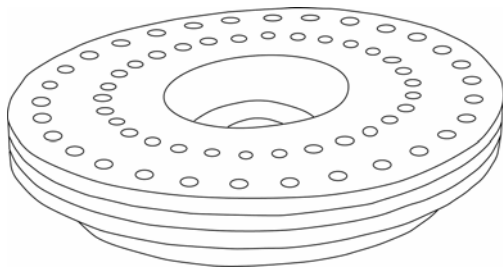

Operation

Select operation mode (“Touch” mode or “On” mode) using the Main Switch

Touch Mode: Unit operates when there is downward pressure on the adapter

On Mode: Unit operates continuously.

Adjust speed using the speed controller. Clockwise to increase the speed.

Turn the main switch to the central “OFF” position when not in use.

Maintenance

Before cleaning, remember to remove the plug from electrical supply.

Periodically, wipe the exterior of the unit using a damp cloth. If necessary, use a non-abrasive and weak detergent solution.

Do not use any chemical cleaners with the unit.

Repair and Servicing

This unit is not designed to be user serviceable.

WARRANTY STATEMENT

LMS guarantees that the device is free from manufacturing defects when shipped from the factory. Should there be any manufacturing defects, LMS will replace the device or repair the defect within 12 months from the date of purchase.


Warranty does not cover units with damaged manufacturer's label, damaged seals, damage through the use of incorrect electrical voltages, and general abuse during usage, transportation or storage. Warranty does not cover parts damaged through normal wear and tear.

The Mixer Uzusio is not designed to be user serviceable. No service manual is available. If the unit is dismantled for any reason, except by the manufacturer or personnel authorized by the manufacturer, all warranties are null and void.

For more information, please contact the distributor below or LMS.

CERTIFICATE OF SALE OF EQUIPMENT

| | |
|-----------------|--|
| Name of company | |
| Address | |
| Model Number | |
| Description | |
| Serial Number | |
| Date of Sale | |

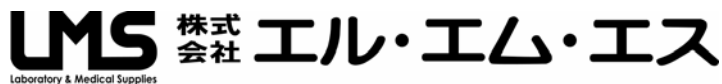

LMS CO., LTD. Tanaka Bldg., 3-6-7, Hongo, Bunkyo-ku,
Tokyo 113-0033, JAPAN

Tel: +813-5842-4171

Fax: +813-5842-4180