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SMALL FUME HOOD GENERAL DESCRIPTION

The small hood comes with a galvanised steel plate case painted with resistant epoxy paint or alternatively with a stainless steel case and it is equipped with a 247 Watt induction motor.

The supporting plane is made of stainless steel.

Under the stainless steel plane there is a service drawer in which tools can be slotted.

The small hood comes with lightening provided by two high performance screw lamps.

The two bag filters supplied with the hood have the following characteristics:

9,200 cm2 nominal filter surface, 20 micron filter nominal capacity.

The control panel comprises the following switches:

- 1 lamp switch 16A
- 1 suction switch 16A
- 1 cleaner switch 16A

...all the switches are covered by a protective rubber shell.

The hood is powered by a 220Volt/50Hz voltage.

The total electrical input is 400 Watt/h (cleaner not included).

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BEHAVIOUR GUIDELINES

1. Use the hood carefully and in a proper way, avoiding any improper or faulty installation breaching this warranty or safety rules in compliance with its manufacturing standards.
2. Concerning hood maintenance, follow very carefully the instructions given in this manual.
3. Before using the appliance, any operator must be perfectly trained on its functions and technical specifications.
4. The hood comes with a one-year period warranty which does not cover any damage or misuse caused by the user.

MAINTENANCE

WARNING! All maintenance must be performed with the unit off and the power plug removed from the outlet. The hood is an easy-to-use appliance requiring little maintenance. Following are a few directions to make it function correctly.

- The bag filters located within the unit must be periodically cleaned, depending on the working hours. However, the cleaning should be carried out any time the sidewalls are saturated with obstructing material and not only when the bag is full.

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PROTECTION GLASS

Always use both hands to adjust the crystal protection glass slant.

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CLEANER SLOT

The cleaner is located on the working top. Above it there is the corresponding power socket.

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LIGHTENING SYSTEM

To replace the lamp, remove the lid on the hood, after having unscrewed the screws. Remove also the second lid.

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After the replacement, reassemble repeating the steps in reverse order.

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POWER SOCKET

The picture on the right shows the cleaner power socket. It also shows the air suction opening.

FILTER MAINTENANCE

To clean or to replace the filter, make sure that the power plug is removed from the socket, then remove the two lids on the hood and extract the bag with the ring nut after having unscrewed the three mounting screws.

The filter can be washed and it must be replaced if the bag is evidently damaged.

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IRIDE INTERNATIONAL S.R.L. Amministratore unico, sig. Enzo Tonin.

DECLARATION OF CONFORMITY

We, IRIDE INTERNATIONAL S.r.l. – furniture for dentist offices and dental laboratories, with registered headquarters in Via San Giuseppe 5 - 33082 Azzano Decimo - Località Tiezzo - PN:

HEREBY DECLARE UNDER OUR OWN RESPONSIBILITY THAT THE PRODUCT “SMALL FUME HOOD” TO WHICH THIS DECLARATION RELATES WERE DESIGNED AND BUILT IN EXCLUSIVE MANUFACTURING IN CONFORMITY WITH THE FOLLOWING HARMONISED STANDARDS:

IEC EN 60601-1(IEC 65-5) of 1991 and its 1994 amendments

IEC EN 60204-1(IEC 44/5 and 44/6)

AND THE REQUIREMENTS OF THE DIRECTIVE:

89/392 of 14/06/1989 and its amendments relating to the harmonisation of the member states' laws for machinery; **89/336/EEC** electromagnetic compatibility; **89/7323/EEC** high and low voltage.

Machinery excluded by Annex 4 of **Directive 89/392/EEC** in compliance with the general requirement **EN 392 CE 89/392 EEC, CE 89/336 EEC, CE 89/7323 EEC, EN292**

IRIDE INTERNATIONAL S.R.L. Governing director, Mr. Enzo Tonin.

IRIDE INTERNATIONAL s.r.l.

Via S. Giuseppe 5 | 33082 Azzano Decimo - PN

tel. +39 0434 633242 | fax +39 0434 632546

welcome@irideinternational.com | www.irideinternational.com