OPERATING MANUAL:  
**PUMP DRIVES**

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A-1299-0818  
Edition 05

## SAFETY TERMS

This manual contains statements regarding safety precautions, which are defined as follows:

**DANGER:** This word and symbol are used in safety messages for hazards that are immediately accessible and capable of causing severe personal injury or death.

**WARNING:** This word and symbol are used in safety messages for hazards that are not immediately accessible, but that present a probability of serious injury if the hazard is not avoided. These safety messages describe precautions that must be taken to avoid the hazard.

**CAUTION:** This word and symbol are used for precautions that must be taken to avoid actions that could damage the equipment.

A **WARNING** or **DANGER** with this symbol indicates probability of serious injury or death exists if an electrical hazard is not avoided.

**Note:** This word is used for information that the operator or technician may use to make the procedure easier to perform.

## SAFETY PRECAUTIONS

**DANGER:** High voltages exist and are accessible in the Cassette Drive. Use extreme caution when servicing internal components.

**WARNING:** Tubing breakage may result in fluid being sprayed from pump. Use appropriate measures to protect operator and equipment.

**WARNING:** Turn drive off before removing or installing tubing. Fingers or loose clothing could get caught in drive mechanism.

## **WARNING: PRODUCT USE LIMITATION**

This product is not designed for, nor intended for use in patient connected applications; including, but not limited to, medical and dental use, and accordingly has not been submitted for FDA approval.

## Introduction

The Manostat CARTER Cassette pump drive provides pump speeds from 3.5 to 200 rpm in 0.1 rpm increments in clockwise or counterclockwise direction.

The PRIME button runs the pump at full speed to prime or purge the tubing.

There are two types of pump operations: Continuous Run Operation and Timer Operation. In Timer Operation, the operator programs the time the pump runs, the time the pump pauses, and the number of cycles in the sequence. Time can be programmed from 0.1 to 59.9 seconds, 1.0 to 999.9 minutes, and 1000 to 9999 minutes.

The digital display shows the pump speed in rpm or percent of maximum speed, counts down timed operation, or shows the number of completed cycles.

The pump may be operated from the control panel (INT) or by remote control by an external current (mA) or voltage (V) supply.

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## Control/Display Functions

Press each button to activate pump function. Press START to run the pump in Continuous Run Operation or Timer Operation. Use up/down (▲, ▼) arrows to change pump speed. Press STOP to stop the pump in Continuous Run Operation or pause in Timer Operation. Press TIME to program automatic-timed settings.

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**A) TIME**—Program automatic running. Sets ON time, OFF time, and number of cycles to be run. Displays number of completed cycles when pump is running in Timer Operation.

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**B) DOWN ARROW (DECREMENT)**—Decreases speed of running pump. Decreases value of a flashing display.

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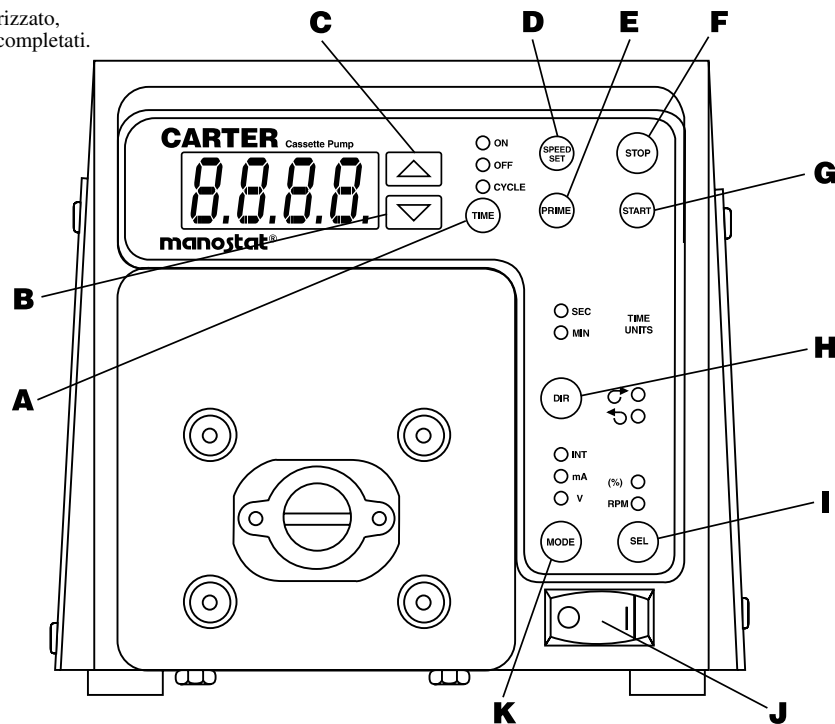
**C) UP ARROW (INCREMENT)**—Increases speed of running pump. Increases value of a flashing display.

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## Setup and Drive Operation

1. Mount cassette pump and load tubing. (See cassette pump manual.)
2. Select motor DIRection (clockwise or counterclockwise).
3. Press SEL to display rpm or % of maximum speed.
4. Select SPEED SET and use ▲ or ▼ arrows to set pump speed.
5. Select MODE (INT, mA, or V).
6. Press and hold PRIME to prime the tubing.

## Purging Tubing

1. Press DIR to select the direction opposite of the pumping action.
2. Press and hold PRIME to purge the tubing.

## Continuous Run Operation

1. Ensure ON, OFF, and CYCLE LEDs are not illuminated. If any are lit, press TIME until all are not lit.
2. Press START to run the pump. Pump runs until STOP is pressed or if MODE is changed.
3. Press ▲ to increase pump speed or ▼ to decrease pump speed.

**NOTE:** Pump restarts automatically after a brownout or power-out condition.

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## Program Timer Operation Function

1. With the pump stopped, press TIME until the ON LED flashes.
2. Use ▲ or ▼ arrows to set the amount of time the pump runs. TIME UNITS displays SEC for seconds or MIN for minutes as the time is being set.
3. Press TIME. The OFF LED flashes.
4. Use ▲ or ▼ arrows to set the amount of time the pump pauses. TIME UNITS displays SEC for seconds or MIN for minutes as the time is being set.
5. Press TIME. The CYCLE LED flashes.
6. Use ▲ or ▼ arrows to set the number of times the pump cycles.
7. Press TIME. The ON, OFF, and CYCLE LEDs illuminate. The pump is ready to operate.
8. Press START to run the pump.

The display counts down the ON time and the OFF time. The ON LED illuminates during the ON time. The OFF LED illuminates during the OFF time. At any time during this operation, press TIME to display the number of completed cycles.

Press ▲ to increase pump speed or ▼ to decrease pump speed.

During Timer Operation, all controls are disabled except for TIME, STOP, and SPEED SET.

Press STOP to pause the operation. Press START to resume the operation.

The speed may be changed at any time during Timer Operation. Press SPEED SET and use ▲ to increase pump speed or ▼ to decrease pump speed.

To cancel the operation, press STOP to pause the pump and then press TIME.

At the end of the programmed number of cycles, the pump stops and remains in Timer Operation.

**NOTE:** The pump does not restart automatically after a brownout or power-out condition. The settings for Time ON, Time OFF, and CYCLE are retained, but the CYCLE count from the last operation is lost. Press START to run the pump.

## Keypad Lockout Enable/Disable

Press and hold SPEED SET. After five (5) seconds, display changes to all dashes. Then, while holding SPEED SET, press PRIME five (5) times.

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## REMOTE CONTROL

Selectable input (0-20 mA, 4-20 mA, 0-10V DC )  
 ±0.5% linearity control  
 2300V isolation potential  
 START/STOP; CW/CCW; PRIME via contact closure (START/STOP only for Timer Operation)

### Remote Control Setup

1. Place the power switch in the off position.

**CAUTION: Power must be turned off before connecting the external remote control cable to prevent damage to the drive.**

2. Connect the cable from the external remote control to the DB15-pin receptacle on the rear panel.
3. Select type of remote control input and output required as follows:
  - a). Press and hold the MODE control while turning the power switch to the on position. After two seconds, release the MODE control. The initial display shows: "inP". After two seconds, the display shows 0-20 or 4-20.

**NOTE:** Press ▲ (increment) or ▼ (decrement) arrows to select 4-20 and 0-20 for current loop control.

- b). Press the MODE control again. The initial display shows: "out". After two seconds, the display shows 0-20, 4-20, or 0-10.

**NOTE:** Press ▲ (increment) or ▼ (decrement) arrows to select 0-20, 4-20, or 0-10 for current loop output.

4. Press the MODE control to select mode of operation. The LEDs indicate the selected mode. Select mA or V.

**NOTES:** During Continuous Run Operation, all functions of the remote control are operable and take precedence over the front panel controls. All contact closures are switch maintained and contact must be held to operate the function. In mA or V mode, the pump does not stop, but runs at the set speed at the completion of a PRIME command.

During Timer Operation, only the START/STOP function is available. Contact closure is momentary. After contact is made and released, the function operates. To stop the function, contact must be made and released a second time.

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## TROUBLESHOOTING AND MAINTENANCE

**WARNING:** Unplug Controller prior to checking or replacing FUSES and/or MOTOR BRUSHES.

### Fuse Replacement

1. Place the power switch in the off position.
2. Disconnect the AC power input line cord from the AC receptacle.
3. Remove and check the fuse and replace if defective.
4. Reconnect the AC power input line cord to the AC receptacle.

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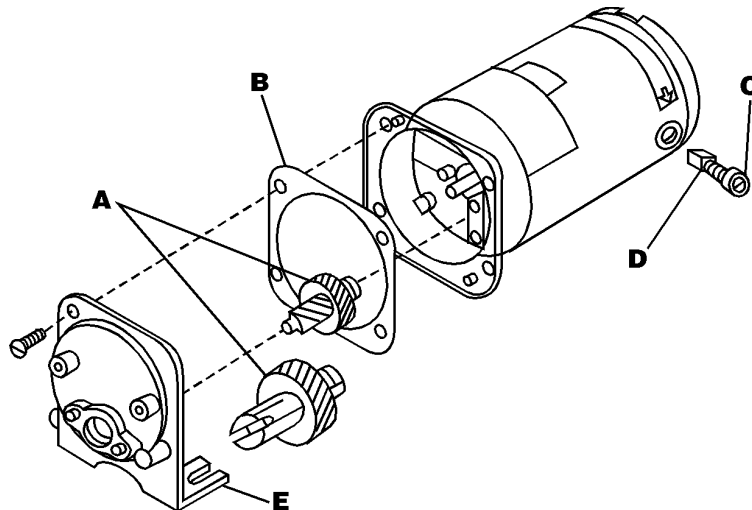
ENGRANAJES

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## MOTOR BRUSH CHECK/REPLACEMENT

**NOTE:** Brushes should be checked every 6 months or 2000 operating hours or if erratic operation occurs.

**DANGER:** High voltages exist and are accessible in the Console Drive. Use extreme caution when servicing internal components.

1. Place the power switch in the off position.
2. Disconnect the AC power input line cord from the AC receptacle.
3. Remove the three screws from each side of the housing and lift off the housing.
4. Carefully unscrew each brush cap on opposite sides of the motor. Withdraw the brush, and examine it for wear.

**NOTE:** One brush is accessible through a hole in the PC board.

**NOTE:** Replace both brushes, if either brush is less than 7.6 mm (0.300 in) long from base to point.

5. Insert brushes and install brush cap on each side of motor.
6. Install housing and secure with the three screws on each side.
7. Connect the AC power input line cord to the AC receptacle.

Contact your dealer if you have service needs.

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### Troubleshooting (cont.)

If an **error message** is displayed, refer to the following list for possible corrective action you can take. If these do not correct the problem, contact your dealer.

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## Cleaning

Keep the drive enclosure clean with mild detergents. Never immerse nor use excessive fluid.

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**REPLACEMENT PARTS  
& ACCESSORIES;  
CLEANING;  
SPECIFICATIONS**

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## SPECIFICATIONS

<b>Output:</b>	
Speed:	3.5 to 200 r/min
Torque:	
Startup	320 oz-in (23 kg•cm)
Continuous	90 oz-in (6.49 kg•cm)
<b>Speed regulation:</b>	
Line	±0.25% F.S.
Load	±0.25% F.S.
Drift	±0.25% F.S.
<b>Display:</b>	Four-digit, seven segment LED
<b>Remote outputs:</b>	Voltage speed output (0–10V DC) Current speed output (0–20 mA or 4–20 mA) Tach output (TTL, 128 to 7680 Hz) Motor running output (N.O. & N.C. contact closure)
<b>Input:</b>	
<b>Supply voltage limits:</b>	
74-000-12131	90 to 130 Vrms @ 50/60 Hz
74-000-12132	200 to 260 Vrms @ 50 Hz
<b>Current, max.:</b>	
74-000-12131	2.3A
74-000-12132	1.2A
<b>Remote Inputs:</b>	START/STOP, CW/CCW, PRIME (Contact closure) Voltage input (0–10V DC) Current input (0–20 mA or 4–20 mA)
<b>Construction:</b>	
<b>Dimensions (L × W × H):</b>	11.2 in × 7.7 in × 7.2 in (284 × 196 × 183 mm)
<b>Weight:</b>	15 lb (6.8 kg)
<b>Enclosure Rating:</b>	IP 23 per IEC 60529
<b>Environment:</b>	
<b>Temperature, Operating:</b>	0° to 40°C (32° to 104°F)
<b>Temperature, Storage:</b>	–45° to 65°C (–49° to 149°F)
<b>Humidity (non-condensing):</b>	10% to 90%
<b>Altitude:</b>	Less than 2000 m
<b>Pollution Degree:</b>	Pollution Degree 2 (Indoor use — lab, office)
<b>Chemical Resistance:</b>	Plastic, gold irridite-coated aluminum, and zinc-plated screws. All materials withstand standard cleaning solvents.
<b>Compliance:</b>	
74-000-12131	UL 508C and CSA C22.2 No. 14
74-000-12132	EU Low Voltage Directive (EN 61010-1) and EU EMC Directive (EN 61326)

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## WARRANTY

Use only MASTERFLEX precision tubing with MASTERFLEX pumps to ensure optimum performance. Use of other tubing may void applicable warranties.

This product is warranted against defects in material or workmanship, and at the option of the manufacturer or distributor, any defective product will be repaired or replaced at no charge, or the purchase price will be refunded to the purchaser, provided that: (a) the warranty claim is made in writing within the period of time specified on this warranty card, (b) proof of purchase by bill of sale or receipted invoice is submitted concurrently with the claim and shows that the product is within the applicable warranty period, and (c) the purchaser complies with procedures for returns set forth in the general terms and conditions contained in the manufacturer's or distributor's most recent catalog.

This warranty shall not apply to: (a) defects or damage resulting from: (i) misuse of the product, (ii) use of the product in other than its normal and customary manner, (iii) accident or neglect, (iv) improper testing, operation, maintenance, service, repair, installation, or storage, (v) unauthorized alteration or modification, or (b) post-expiration dated materials.

THIS WARRANTY IS THE EXCLUSIVE REMEDY OF THE PURCHASER, AND THE MANUFACTURER AND DISTRIBUTOR DISCLAIM ALL OTHER WARRANTIES, WHETHER EXPRESS, IMPLIED, OR STATUTORY, INCLUDING WITHOUT LIMITATION, WARRANTIES OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE. NO EMPLOYEE, AGENT, OR REPRESENTATIVE OF THE MANUFACTURER OR DISTRIBUTOR IS AUTHORIZED TO BIND THE MANUFACTURER OR DISTRIBUTOR TO ANY OTHER WARRANTY. IN NO EVENT SHALL THE MANUFACTURER OR DISTRIBUTOR BE LIABLE FOR INCIDENTAL, INDIRECT, SPECIAL OR CONSEQUENTIAL DAMAGES.

*The warranty period for this product is two (2) years from date of purchase.*

## PRODUCT RETURN

To limit charges and delays, contact the seller or manufacturer for authorization and shipping instructions before returning the product, either within or outside of the warranty period. When returning the product, please state the reason for the return. For your protection, pack the product carefully and insure it against possible damage or loss. Any damages resulting from improper packaging are your responsibility.

## TECHNICAL ASSISTANCE

If you have any questions about the use of this product, contact the manufacturer or authorized seller.

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**GARANTIE  
GARANTIE  
GARANTIA  
GARANZIA**

NOTES

### Operating Manual Addendum Cassette Loading Capacity

Use the chart below for maximum cassette loading.

Soft material is silicone. Stiff materials: PHARMED®, SANTOPRENE®, TYGON®, PVC, and VITON® fluoroelastomer. N/R = not recommended. CARTER's continuous output torque is 90 oz-in. (6.5 kg•cm).

Tubing Link (mm ID)	Tubing Material	Max # of Cassettes	Tubing Link (mm ID)	Tubing Material	Max # of Cassettes
<b>74-048-00000</b>			<b>74-046-00000</b>		
.19	Soft	4	13	Soft	2
	Stiff	4		Stiff	2
.25	Soft	4	14	Soft	2
	Stiff	4		Stiff	2
.89	Soft	4	16	Soft	2
	Stiff	4		Stiff	2
1.42	Soft	4	25	Soft	2
	Stiff	4		Stiff	1
2.06	Soft	4	17	Soft	2
	Stiff	4		Stiff	N/R
2.79	Soft	4	18	Soft	2
	Stiff	4			
<b>74-128-00000</b>			<b>74-126-00000</b>		
.19	Soft	12	13	Soft	6
	Stiff	12		Stiff	6
.25	Soft	12	14	Soft	6
	Stiff	12		Stiff	6
.89	Soft	12	16	Soft	6
	Stiff	12		Stiff	6
1.42	Soft	12	25	Soft	6
	Stiff	12		Stiff	4
2.06	Soft	12	17	Soft	4
	Stiff	10		Stiff	3
2.79	Soft	12	18	Soft	3
	Stiff	7			
<b>74-083-00000</b>			<b>74-083-00000</b>		
.19	Soft	N/R	13	Soft	4
	Stiff	N/R		Stiff	4
.25	Soft	N/R	14	Soft	4
	Stiff	N/R		Stiff	4
.89	Soft	8	16	Soft	4
	Stiff	8		Stiff	3
1.42	Soft	8	25	Soft	4
	Stiff	8		Stiff	2
2.06	Soft	8	17	Soft	4
	Stiff	8		Stiff	2
2.79	Soft	8	18	Soft	4
	Stiff	8			
<b>74-084-00000</b>			<b>74-084-00000</b>		
.19	Soft	N/R	13	Soft	4
	Stiff	N/R		Stiff	4
.25	Soft	N/R	14	Soft	4
	Stiff	N/R		Stiff	4
.89	Soft	8	16	Soft	4
	Stiff	8		Stiff	3
1.42	Soft	8	25	Soft	4
	Stiff	8		Stiff	2
2.06	Soft	8	17	Soft	4
	Stiff	8		Stiff	2
2.79	Soft	8	18	Soft	4
	Stiff	8			

These ratings are a guideline only. Please consult the factory if you need to exceed these ratings.

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