technotrate

Manuale operativo

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Operating Manual

Please study carefully and follow step by step !

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Safety Instructions

1. Every user must study these instructions carefully before using the instrument.
2. General operation and safety instructions must be followed. Eye protection, protective clothing and glove must be worn.
3. The instrument must only be used for intended purpose and utilized within the component tolerances of material resistance.
4. Always work with the instrument so as not to endanger the operator or other persons.
5. Do not apply force to the instrument.
6. Do not drop the instrument.
7. Do not immerse the instrument in liquids.
8. Only use original spares and accessories.
9. Repairs must only be carried out by trained personnel. Do not open the instrument. Such actions will nullify the guarantee!
10. Do not use the instrument in explosive environments.
11. Do not expose the instrument to temperatures exceeding 50°C. Do not autoclave !
12. Keep the instrument away from fire or flames.

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Index

Safety Instructions	2
Use	4
Commissioning of the instrument	4
Installation of Discharge tube	5
Installation of Suction tube	5
Seating onto the flask	5
Charging	6
Switch on	6
Using the burette	6
Purging	6
Suction	7
Titration	8
Dosing	9
Draining	10
Storage of data	10
Calibration mode	12
Common instructions to user	13
Cleaning	13

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Use

The technorate Digital Burette is intended to be used for titration and dosing within following limits:

15 - 40°C of reagent and instrument

When using according to the instructions, only these parts of the Digital Burette will come into direct contact with following chemically resistant material :

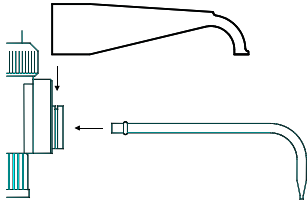
- FEP (Suction hose, Discharge tube)
- Tefzel® (Valve block)
- Borosilicate glass (Valves, KPG Cylinder)
- Ceramic, Platinum-Iridium (Valves)

Never use the instrument with

- Hydrofluoric acid
- liquids which corrode PTFE, FEP, Tefzel®, Platinum-Iridium or where catalytic changes to these material can occur
- crystallizing, particulating or precipitating solutions (e.g. suspensions)

the adapter included with the supply are of polypropylene and may only be used with non PP-aggressive media!

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
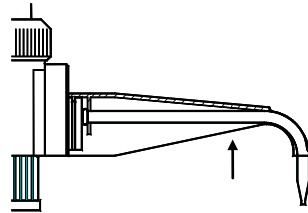
Installation of Discharge tube

2

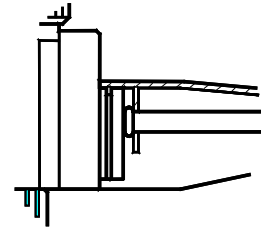Put supplied Discharge tube into the existing opening and clip-on the protective sleeve (1). Fix the Discharge tube by pressing slightly into the protective sleeve (2). The Discharge tube must be installed that way so that the bulge is placed between valve body and fixing bridge of the protective sleeve (3).

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Installation of Suction tube

Cut suction tube and reverse dosing tube to desired length and put into the sockets of the bottom part of the instrument. Check for firm seating.

Seating onto the flask

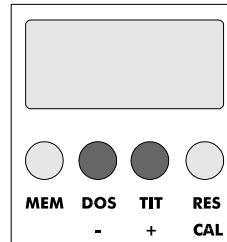
Seat instrument onto the flask, turn to desired position to flask and screw on tightly by simultaneous seizing of the burette and turning of the GL45 cap.

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Charging

If necessary, charge the built-in NiCd batteries with the supplied charging unit. First put the charging unit into mains socket and then the plug into the bush at the back of the burette.

Switch on

Switch on the instrument by pressing one of the buttons *DOS* and *TIT*. While pressing it appears „dos“ or „tit“ on the display.

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

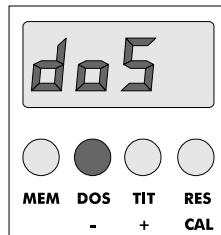
2

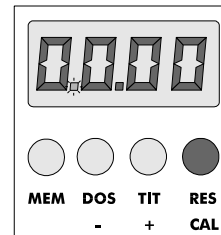3

Turn selector switch to position *discharge* (1) and press button *DOS* (2). While pressing the button *RES* adjust the display in the dosing mode to 00.00 (3). Turn handwheels in an even but not too fast speed upwards until the required volume appears in the display.

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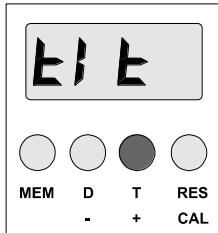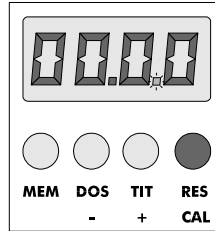
Titration


2

Operate TIT switch (1) (in the display flashes the dot between the third and fourth digit) and by pressing the button RES (2) adjust the display in the titration mode to 00.00. Place a test flask under the discharge tube and turn the handwheels downwards without interruption and not exceeding the titration speed of 1.5 ml/sec.


The titrated volume can be read off immediately after reaching the tit point (e.g. colour change) (3).

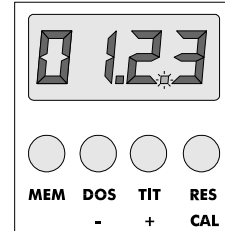For further titration simply adjust display to 00.00 by pressing the button RES and proceed with titrating. In case the burette cylinder is empty before the titration is finished, turn handwheels upwards again in order to suck additional titration liquid out of the flask. If necessary control the amount of liquid by pressing the buttons DOS (the dot between the first and second digit of the display flashes) and RES. The display of titrating mode is not influenced. Carry on with titration after pressing the button TIT.

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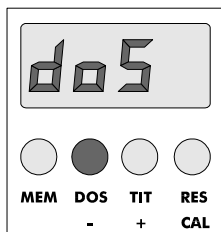
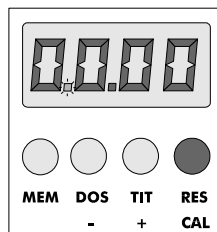
Dosing

2

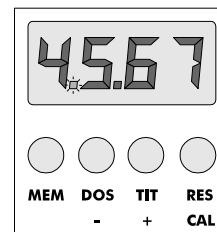
The technorate digital burette can be also used for dosing of suitable media. Press the button *DOS* (1) (in the display flashes the dot between the first and second digit), adjust the display to 00.00 by pressing the button *RES* (2) and suck the amount of medium to be dosed (3). Place a suitable receptacle under the discharge tube and then discharge the entire amount by turning the handwheels downwards.

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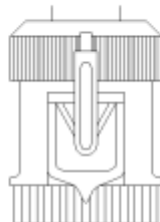Draining

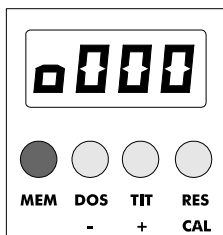
To drain the discharge tube without losses turn the selector switch to position *draining*. The remaining liquid in the discharge tube flows back into the flask. This position of the selector switch does not allow titrating or dosing by mistake because the discharge of liquids is prevented. If liquid erroneously has been sucked from the flask, turn selector switch to position *draining* and dose back liquid into the flask.

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
Storage of data


The technorate possesses the facility to store in the chip up to 500 values which after the completion of e.g. serial titration can be transferred by WITEGSTAT software using WITOFACE modem and processed on any conventional PC. For details refer to the instruction manual supplied with WITOFACE.

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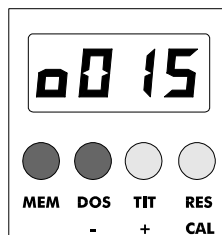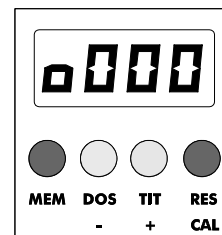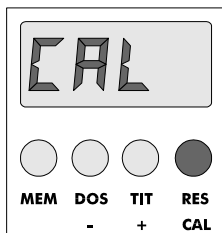
For the storage of measured values, press button *MEM* (1). It appears while the button is depressed the display „o000“ indicating the first stored value. Every further press on the button *MEM* will store a value and the display of the stored values will increase by 1.

If the *MEM*-button kept depressed and simultaneously pressing one of the buttons *DOS* or *TIT* (2), the data storage decreases by 1 for every press, this means that the last value each is erased. To erase the total storage, depress *MEM*-button and simultaneously press button *RES* (3). It appears o000 in the display. The storage is empty and can again store up to 500 values.

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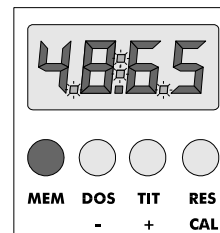
Calibration Mode

The instrument is factory calibrated to 20°C aqueous solutions. For solutions of varying density and viscosity, the technotrate can be adjusted by the user in case of changed conditions provided a laboratory balance with a resolution of <1 mg is available.

- 5 titrations of nominal volume of 50.00 ml to be weighted out on the balance and applying the temperature correction factor (our testing instructions PA...) the true values to be investigated
- Average the 5 resultant values
- Depress buttons RES/CAL until in the display appears „CAL“ (1), release button and in the display the dots flash between the first and second as well as the third and fourth digit.
- Adjust the display (e.g. 50.00) with the buttons DOS/- and TIT/+ to the final average value (2)
- To activate the calibration value press button MEM (3) (in the display a colon flashes additionally between the 2nd and 3rd digit)
- Press button „RES/CAL“ to return to operation mode

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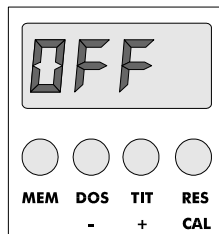
- During the work with this calibration value, in the display a colon flashes between the second and third digit
- In order to return to the factory settings, depress button „RES/CAL“ until in the display „CAL“ appears, then deactivate the calibration value by pressing the button *MEM* and return to operation mode via pressing „RES/CAL“. The flashing of the colon on the display disappear

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Common instructions to user

In case the burette remained for longer period screwed to the flask and not in regular use, we recommend to purge the instrument prior to again titration or dosing.

Caution!! The instrument is not equipped with an „off“ switch“.

The technorate automatically is switching off after 3 minutes of deactivation. The display of the instrument indicates „OFF“. To reactivate simply press button „DOS“ or button „TIT“. In the display appears now each the last titration and dosing values.

Cleaning

The instrument should be cleaned regularly but always on

- reagent change
- long storage
- use of crystallizing media

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Warning !

When cleaning observe all safety instructions. Components of the instrument are contaminated with titrating liquid! Suitable protective clothing, safety glasses and gloves must be worn!

- Drain instrument and place it along with the bottle inside a washbasin
- Unscrew instrument from the bottle and pull out until suction and reverse dosing tubes are no longer immersed
- Turn selector switch to position *purging* and move handwheels several times upwards and down wards until the suction and reverse dosing tubes are free from liquid
- Immerse tubes in suitable cleaning solvent and rins thoroughly by applying suction and reverse dosing; turn selector switch to *discharge* and repeat operation in order to clean discharge tube and discharge channels

Finally rinse with distilled water

Accessori	Accessories	Specifiche	Art.-n.
Interfaccia Witolink Kit die collegamento technorate-PC compressa interfaccia, software e fili	Witolink technorate-PC-Connection Kit incl. interface, software, wires		83022
Bottiglia	bottle	1000 ml, GL45	83011
Bottiglia	bottle	2500 ml, GL45	83012
Cannula di scarico technorate	technorate discharge tube		83019
Tubo aspirazione technorate	technorate suction tube	300 mm	83015
Tubo di circolazione	Circulation tube	300 mm	83020
Carica-batteria technorate	technorate battery charger	220 V	83021
Set di 5 adattatori	Set of 5 adapters	A25, 28, 38, 40, 45	83023
Tubo cloruro di calcio	Calcium Chloride tube		83024
Valvola di controllo	Check valve		83026