
**Operating instructions  
Bedienungsanleitung  
Mode d'emploi  
Instrucciones de manejo  
Istruzioni d'uso**

**METTLER TOLEDO  
LC-P43 Printer  
Drucker LC-P43  
Imprimante LC-P43  
Impresora LC-P43  
Stampante LC-P43**

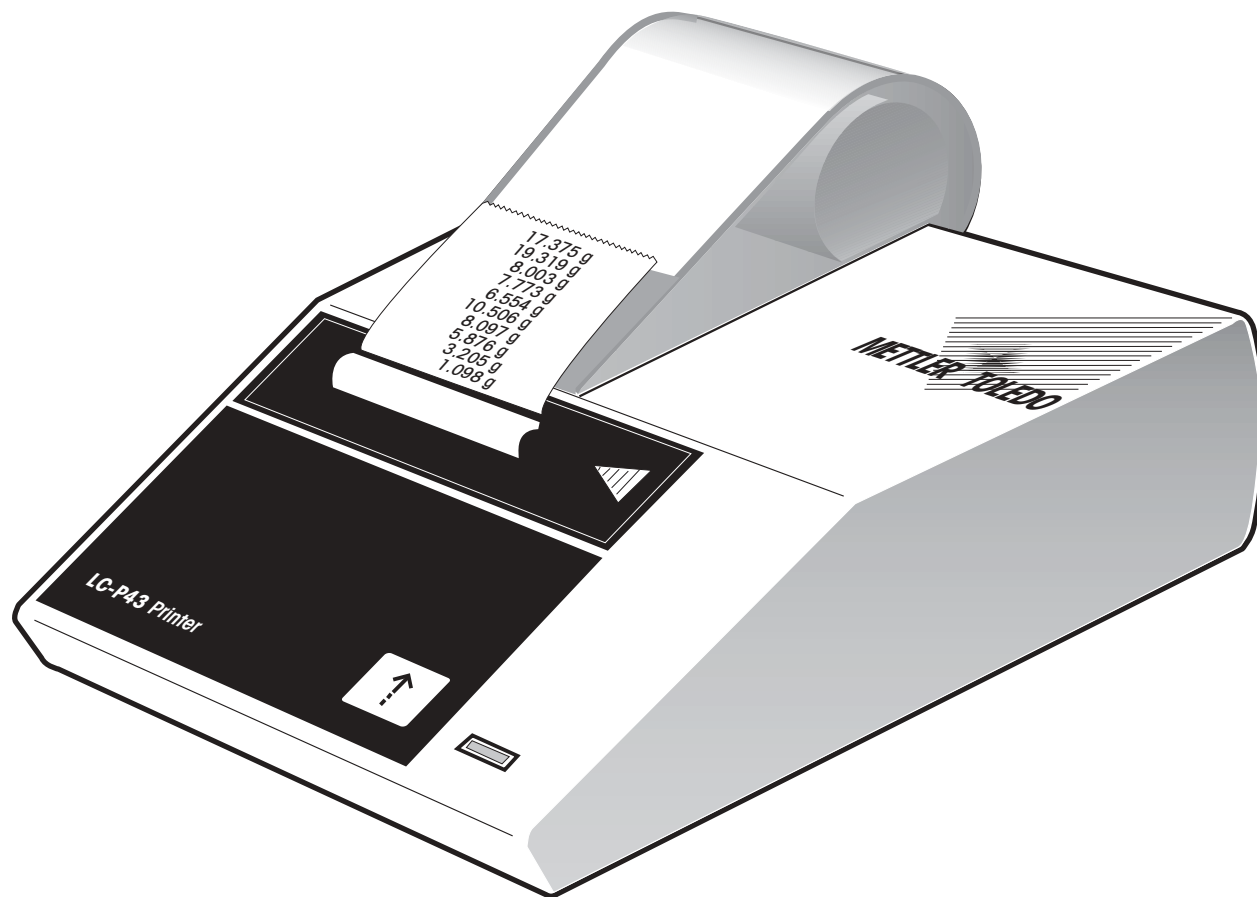**METTLER TOLEDO**

## Overview, Übersicht, Sommaire, Indice, Indice

Introduction .....	4	Einführung .....	11	Introduction .....	18	Introducción .....	25	Introduzione .....	32
Startup procedure .....	4	Inbetriebnahme .....	11	Mise en service .....	18	Puesta en funcionamiento .....	25	Messa in servizio .....	32
Operation .....	5	Bedienung .....	12	Utilisation .....	19	Manejo .....	26	Impiego .....	33
Printer test .....	6	Test .....	13	Test .....	20	Test de impresora .....	27	Test della stampante .....	34
Maintenance .....	7	Wartung .....	14	Entretien .....	21	Mantenimiento .....	28	Manutenzione .....	35
Malfunctions .....	7	Störung .....	14	Pannes .....	21	Fallos .....	28	Disturbi .....	35
Maintenance (Service) .....	8	Wartung (Service) .....	15	Maintenance (Service) .....	22	Mantenimiento .....	29	Manutenzione .....	36
Cleaning .....	8	Reinigung .....	15	Nettoyage .....	22	Limpieza .....	29	Pulizia .....	36
Accessories, Spare parts ...	9	Zubehör, Ersatzteile .....	16	Accessoires, Pièces de rechange .....	23	Accesorios, repuestos ...	30	Accessori, ricambi .....	37
Technical data .....	9	Technische Daten .....	16	Caractéristiques techniques .....	23	Características técnicas ..	30	Caratteristiche tecniche ...	37

**English Page 4**

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**Español Page 25**

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## Introduction

The LC-P43 is a rugged and simple to operate dot matrix printer for balances from METTLER TOLEDO with a LocalCAN universal interface (balance lines AB, PB, CB, GB, College-B, SB with LC-B option, AG, PG, SG, PR, SR and PG-S with LC-G option). The printout on normal paper fulfills the requirements of modern quality assurance systems (GLP, GMP, ISO 9000, etc.). These define, among other things, the principles and duties for the user of measuring instruments or inspection, measuring and test equipment. The LC-P43 complies with the principle of traceability of measurements by precisely documenting the origin of data, as well as calibration procedures. It fulfills the obligation to archive measurement results by its chemically resistant, thermally stable and lightfast print. The LC-P43 prints all data sent by the balance without restriction, that is not only weight values and calibration procedures, but also naturally the results of all applications such as piece counting built into the balance.


## Startup procedure

### Cautionary notes

For reasons of safety, the housing of the printer must not be opened by the user. If technical faults appear or if the printing unit or power line fuse has to be changed, please notify METTLER TOLEDO service.


- The LC-P43 Printer may be used only indoors and on no account in hazardous areas.
- The electronics of the LC-P43 Printer are protected against the ingress of splashing water and dust. However, as the printer housing is not tight the printer should not be used in the presence of liquids.
- Comply with cleaning instructions (see page 8), protect power plug against wet conditions.
- Never open the printer housing. Should you ever have problems with your printer, please contact the responsible METTLER TOLEDO dealer.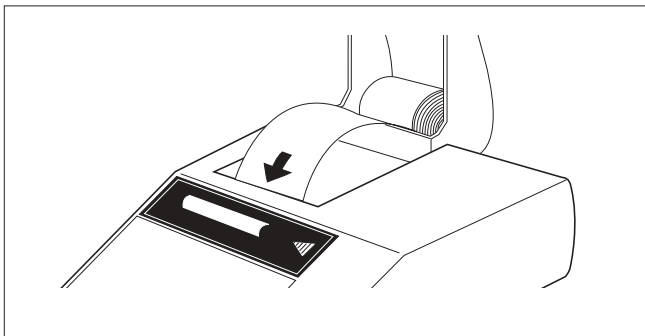


### Inserting paper


- Feed paper through the slot in the printing unit in the direction of the arrow.
- Switch on printer with toggle switch at the rear.
- Press and hold  until sufficient paper has been fed in

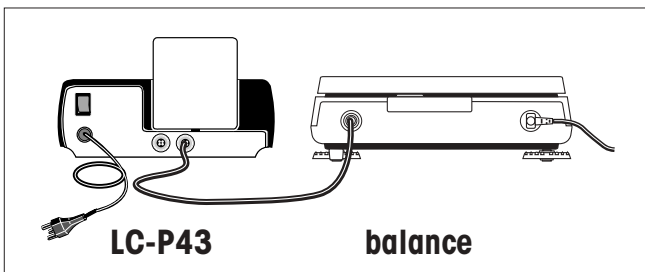### Attaching printer, printing

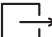- Attach data cable of the LC-P43 to the data interface of the balance.  
The printer is now ready for operation!
- Pressing the balance key with the  symbol prints out the weight value.  
(PR- and SR-balances: weight with time, date and sample number.)

## Operation

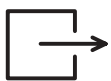
- Pressing the balance key with the  symbol prints out the stable weight value.  
As long as the balance does not show a stable weight value, nothing will be printed out.  
(With PR- and SR-balances, an automatic printout is also possible.)

- Pressing the printer key  advances the paper 1 line (line feed).

## Printer test

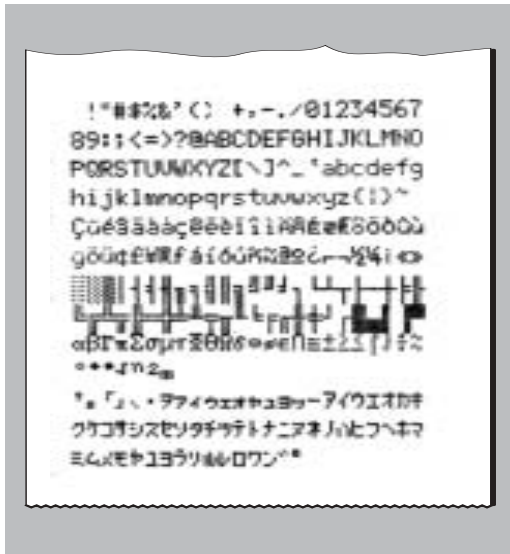The LC-P43 is equipped with a self-test which automatically prints out the character set of the printer.


### Start the test

- Switch on LC-P43 and at the same time press and hold  for a few seconds.

### Terminate the test

- Switch off the LC-P43.

## Maintenance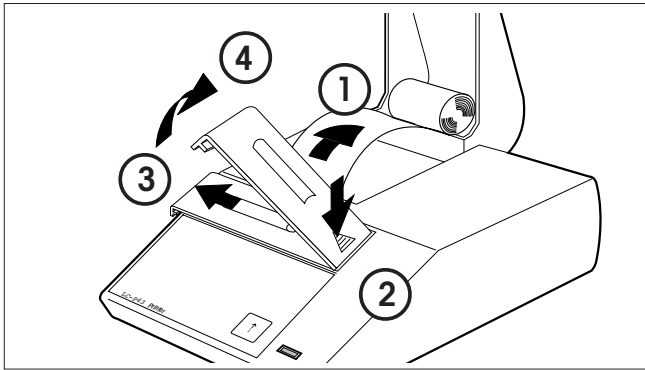


### Changing paper

- Insertion of a new roll of paper is described on page 5. Order number for paper, see page 9.

### Replacing the ribbon

- 1 Pull paper out of printing unit.
- 2..4 Remove ribbon cover.
- 5 Remove ribbon cartridge (press in direction of arrow).
- Insert new ribbon (if need be, take up slack with tensioning wheel 6).
- Refit ribbon cover.
- Insert paper and feed through slit in the printing unit as far as it will go.
- Press and hold  until sufficient paper has been fed in.

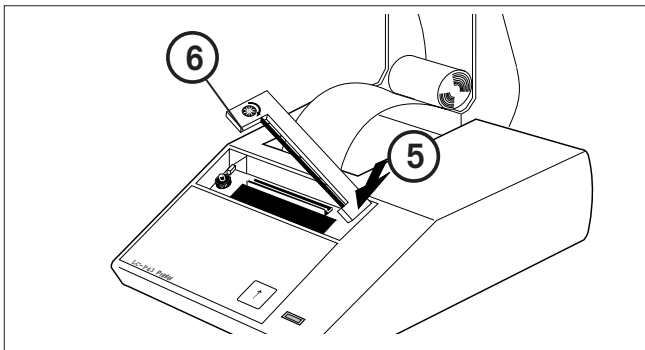## Malfunctions

### Malfunction/Message

Printout unreadable

Green pilot lamp does not light up  
Green pilot lamp flashes slowly

### Possible cause

- Ribbon worn/entangled
- Service life of printing unit at an end
- No power, fuse defective
- Balance waiting for stability/cable not attached

### Rectification

- Change/tension ribbon
- Please contact METTLER TOLEDO service
- Please contact METTLER TOLEDO service

## Maintenance

### Service

Under normal ambient conditions, the LC-P43 Printer requires practically no maintenance thanks to its rugged construction. The METTLER TOLEDO service department is available should servicing be required. Please ask your METTLER TOLEDO dealer for details.

## Cleaning

As the printer housing is made of top grade, resistant materials, all commercially available cleaning agents may be used.

<b>Accessories, Spare parts, expendables</b>
--

Accessories, spare parts, expendables	Type/Order number	Standard equipment
Paper roll, set of 5	72456	2
Ribbon cartridge, black, minimum order quantity 2	65975	1
Extension cable/connection cable LocalCAN 2 m	LC-LC2 cable	–
Extension cable/connection cable LocalCAN 5 m	LC-LC5 cable	–
T-socket for devices with only one LocalCAN socket	LC-LCT	–
Spare printing unit (may be changed only by METTLER TOLEDO service)	–	–
Power line fuse 230 V (slow-blow 100 mA, 250 V) + 115 V (slow-blow 200 mA, 250 V) (may be changed only by METTLER TOLEDO service)	–	–

<b>Technical data</b>
-----------------------

<b>Printer</b>	
Printing unit	dot matrix printer 5x9 dots, 24-character line length, Epson IBM character table No. 4
Printing speed	min. 1 line per second
Ribbon cartridge	exchangeable, black
Paper roll	normal paper, 58 x dia. 51 mm, integrated in housing, commercial size
<b>Printer functions</b>	Printing, lightfast and thermally stable (GLP, GMP, ISO9000), line feed, self-test.
<b>Interface</b>	<b>LocalCAN</b> , general-purpose interface with 2 equal priority outputs for the simultaneous attachment of up to five different peripherals (e.g. auxiliary display, foot switch, printer; computer via RS-232-C).

**Designed test conditions**

Power line voltage, frequency	115 V or 230 V, admissible fluctuations -20%+15%, 50/60 Hz
Power consumption	140 mA or 70 mA
Altitude	4000 m
Temperature range	0...40 °C
Humidity	max. relative humidity 80% for temperatures to max. 31 °C linear decreasing to 50% at 40 °C.
Installation category	II
Pollution degree	2
Dimensions/weight	W x D x H = 157 x 210 x 85 mm / net 1.5 kg (incl. paper roll)

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
**To protect your METTLER TOLEDO product's future:  
METTLER TOLEDO Service assures the quality, measuring accuracy and preservation  
of value of all METTLER TOLEDO products for years to come.  
Please send for full details about our attractive terms of service.  
Thank you.**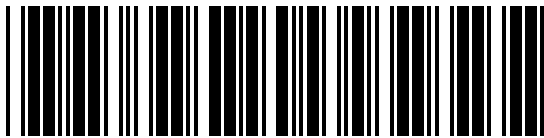

\*P705177\*

Subject to technical changes and to the availability  
of the accessories supplied with the instruments.

Printed on 100 % chlorine-free Paper,  
for the sake of our environment.

© Mettler-Toledo GmbH 1998      705177C Printed in Switzerland 9806/2.45

**Mettler-Toledo GmbH, Laboratory & Weighing Technologies**, CH-8606 Greifensee, Switzerland  
Phone +41-1-944 22 11, Fax +41-1-944 30 60, Internet: <http://www.mt.com>