StarFireST

Operator's Manual
Mode d'emploi
Istruzioni per l'uso
Bedienungsanleitung
Manual del Operador

ARGOS
TECHNOLOGIES *manufacturing solutions for science*

Argos Technologies, Inc.
1551 South Scottsdale Court
Suite 200
Elgin, IL 60123
Toll Free: 800-886-8675
International: 847-622-0456
www.argos-tech.com

Table of Contents

Contents	2
Introduction	3
Hazard Warnings & Caution Notes	3
Hazard Warnings used in this instruction manual	3
Hazard Warnings used on the Device	3
Important Safety Information	4
Safety Features of StarFire ST	4
Using StarFire ST	4
General Safety Notes	5
Restrictions for use	5
Warranty Information	6
Repair Service	6
Product Description	7
Package contents	7
Front View	7
Rear View	7
Burner Head	8
Bottom View	8
Getting Started	9
Connecting StarFire ST to a gas supply	9
Changing the nozzles	9
Fitting the appropriate gas nozzle	9
Installing and removing adapters for gas connection	10
Connecting the gas cartridge	11
Removing and changing the gas cartridge	11
Setting the right flammable mixture	12
Connecting the foot switch	12
Removing and attaching the burner head	12
Connecting StarFire ST to a power supply	12
Operating StarFire ST	13
Igniting the flame	13
Using the ignite button	13
Using the footswitch	13
Working with liquids	13
Maintenance, Care and Safety Precautions	14
Care and Maintenance	14
Troubleshooting (FAQ)	15
Technical Data	16
Accessories and Spare Parts	17

Introduction

Thank you for your recent purchase of the Argos StarFire ST, the safe alternative to the standard Bunsen Burner.

This instruction manual contains all the information required to safely use and maintain the StarFire ST.

Hazard Warnings & Caution Notes

Hazard Warnings used in this instruction manual

**DANGER**

Indicates possible danger of personal injury or death. It is important that you comply with the instructions and precautions.

**CAUTION**

Indicates possible damage to the instrument. Comply with the instructions.

**NOTE**

Provides important notes on the correct operation of StarFire ST.

Hazard Warnings used on the Device

**ATTENTION HOT**

Risk of burns from the flame.

Important Safety Information

StarFire ST is an automatic Bunsen burner that can be used to heat or sterilize samples, flasks or other laboratory materials. Unlike traditional Bunsen burners, StarFire ST is not intended for continuous heating applications.

Always use this product in accordance with these instructions. Where appropriate, safety goggles and protective clothing may be required for your safety.

Safety Features of StarFire ST

- If the flame is not ignited within 5 seconds, the flame detection will shut off the gas supply.
- If the temperature exceeds the safe operating level the gas supply to the unit is automatically shut off.

Using StarFire ST

StarFire ST has been designed to incorporate several safety features that comply with industry recognized safety standards. The unit should only be operated by trained personnel following the operating instructions detailed in this manual.

StarFire ST should never be used in disrepair or operated by untrained personnel. Doing so may result in potential safety risks to the laboratory.

To ensure the safe operation of the StarFire ST, all personnel that operate the burner must have read and understood the manual. Should you have any questions regarding its operation please contact our Technical Support department prior to operation.

Contact your local safety officer if you have any questions about local or corporate regulations and guidelines related to the safe operation of our instrument.

Important Safety Information (cont.)

General Safety Notes

- Never leave an activated StarFire ST unattended
- When unpacking, be sure to check for any transportation damage. Do not operate if damage is detected.
- Always remove the gas cartridge *before* removing the gas cartridge adapter.
- There are no serviceable parts inside the StarFire ST, contact Argos Technologies or a trained service dealer should your instrument need repair or service.
- Do not change or alter the device in any manner. Doing so nullifies the warranty and may pose a safety risk.
- Only use Argos parts, supplies and accessories with the StarFire ST unit.
- Always wear proper safety equipment (especially protective eyewear) when operating the StarFire ST.
- StarFire ST should only be operated in well ventilated areas and away from any flammable or combustible materials.
- To ensure safe operation, all gas connections should be tightened and checked for leaking.
- Only approved safety gas tubing should be used with the StarFire ST.
- Keep hands and face away from the burner head at all times.
- The burner head remains hot after the flame is extinguished. Do not touch. Allow sufficient time for the head to cool down before transporting, servicing or cleaning the unit.
- Do not pour or spill materials into the burner head. Doing so may prevent the unit from operating properly.

Important: You must turn off the gas supply when:

- The unit is initially installed or moved to a new location
- The gas cartridge is replaced
- During extended periods when the unit is not in use
- Any time you clean or service the unit

Restrictions for use

CAUTION

The StarFire ST is not intended to be operated by turning the unit upside down. Never carry the StarFire ST by the burner head as the unit may detach.

Warranty Information

Your Argos StarFire ST is warranted to be free of defects in materials and workmanship for a period of 12 months from the date of purchase, providing that:

- The unit has not been opened or attempted to be opened by the end user
- The user ensures that it has been carefully cleaned and decontaminated
- The unit has been operated according to these instructions and not in any improper way

In order to qualify for warranty repair; the customer must return the warranty card supplied with the unit that includes all of the information required to validate proof of purchase. Should warranty service be required, repair or replacement will be provided by Argos without charge. Freight shipping the unit to Argos will be the responsibility of the purchaser. There are no other warranties of application or merchantability.

Repair Service

Should your StarFire ST require service or repair, please contact Argos Technologies Customer Service. They will issue a Return Authorization Number and advice on the proper packing and method of shipment.

Customer Service:

Within US: 800-886-8675

Outside US: 847-622-0456

Argos Technologies, Inc.
1551 South Scottsdale Court
Suite 200
Elgin, IL 60123
Attn: Technical Services

Be sure that your RA number is on the shipping label.

Please use the same above numbers if you require any technical assistance or support.

Product Description

Package contents

- StarFire ST Burner
- Universal Power Supply
- UK Adapter
- US Adapter
- EU Adapter
- Alternate Nozzle
- Warranty Registration Card
- Instruction Manual
- 10mm gas adapter for mains gas supply
- Footswitch

Front View

1. Burner Head
2. GAS KNOB: to adjust the gas supply (flame size)
3. AIR KNOB: to adjust the air supply (flame temperature)
4. IGNITE BUTTON: to switch the device on and off. Ignites flame


Rear View

5. Foot Switch connection (jack)
6. Mains plug socket (mains plug)
7. Gas adapter opening (for engaging)


Product Description (cont.)

Burner Head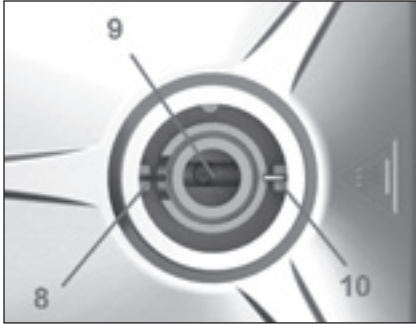

- 8. Ignition Electrode
- 9. Burner chamber with flame opening
- 10. Electrode of the flame monitor

Rear View

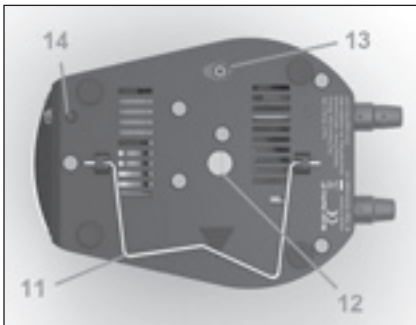- 11. Folding Stand, to position device at an angle
- 12. Brass screw, access to the nozzle
- 13. Alternative Nozzle
- 14. Release button for adapters for gas connection

Getting Started

Please ensure that all packaging materials have been removed before you start using your StarFire ST. Place StarFire ST on a flat and dry surface.

Connecting StarFire ST to a gas supply

StarFire ST can be used with natural gas (gas mixture, >85% methane, further alkanes) and commercial butane/propane gas (gas mixture, 65%-10% butane, 20%-40% propane). When connecting to natural gas be sure to use only approved and tested gas safety hose with an inner diameter of 10 mm.

Changing the nozzles

To exchange gas nozzles, follow the following procedures:

- Stand the unit upright to access the bottom of the case.
- Remove the brass screw (1) from the case bottom using a flat head screwdriver
- Tilt or turn the device over until the nozzle drops out of the burner head assembly.
- Locate the nozzle holder next to the brass screw that holds the extra nozzle.
- Remove the nozzle from the tube and insert the new nozzle into the burner head assembly.
- Replace the brass screw and tight to the stop point.

Fitting the appropriate gas nozzle

The E nozzle for natural gas is fitted ex-factory for StarFire ST. Should you plan to use either propane or butane as fuel, please change the nozzle.

Type of Gas	Gas Input Pressure	Inscription on Nozzle
Natural gas (methane)	20 mbar	E
Butane/propane gas	50 mbar	P

CAUTION

StarFire ST must only be operated with the correct nozzle that corresponds to the type of gas used. Only use nozzles supplied with StarFire ST. While changing nozzles the user should inspect the o-ring located in the brass screw (see above) for damage and replace if necessary.

Getting Started (cont.)

Installing and removing adapters for gas connection


StarFire ST is supplied with both a barbed 10mm tubing adapter and a ¼" left handed threaded adapter. Choose the adapter that is appropriate for your laboratory setting. These adapters are to be used when using house or natural gas and not intended for use with the CV360 gas cartridge.

- Insert the appropriate adapter in the gas adapter opening of StarFire ST, until it clicks into position.

CAUTION

When changing the gas adapter ensure the device is switched off and disconnected from the mains supply, and the gas supply is shut off.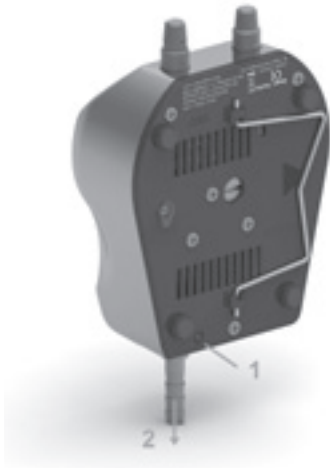

- To remove the adapter, use a pointed tool, such as a ball point pen, to press the adapter release button on the underside (1) of the device and pull the adapter away from the device (2)
- Insert the new adapter into the gas adapter opening until it clicks into position.

Getting Started (cont.)


Connecting the gas cartridge

WARNING

To operate the StarFire ST using a gas cartridge, the P nozzle must be used in the burner head. Only use gas cartridges between ambient temperatures of 15-35°C.

The CV 360 cartridge is connected as follows: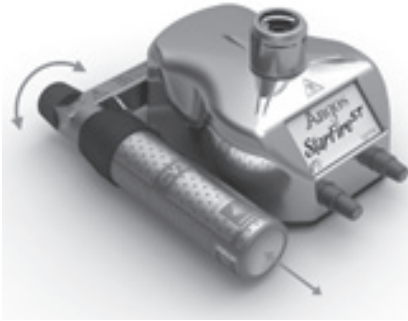

- Insert the gas cartridge adapter in the adapter opening until you hear a “click”
- Insert the gas cartridge into the cap until it audibly clicks into position
- The gas cartridge is now inserted
- Turn the black knob fully in the “ON” direction to allow the gas to flow.
- The flame can be ignited using any of the two operating modes with the gas canister connected.
- Verify that there are no leaks prior to engaging flame.

CAUTION

Care must be taken not to dislodge the gas canister during use.

Removing and changing the gas cartridge

WARNING

Do not remove or replace the gas cartridge unless the gas adapter is connected to the device and shut off (by turning the black knob in the “OFF” direction till it stops)

The CV 360 gas cartridge is removed and changed as follows:

- The gas adapter should be attached to StarFire ST
- Turn the black knob in the “OFF” direction until the gas is shut off
- Pull the gas cartridge away from the gas adapter (towards the front of the device)
- Push a new gas cartridge into place (see “connecting the gas cartridge” above)
- When pulling the cartridge from the adapter small amounts of gas may leak out but are not considered dangerous

Getting Started (cont.)

Setting the right flammable mixture

You can adjust the size of the flame by rotating the GAS KNOB on the front of the unit. Rotating the knob reduces the gas supply and lowers the flame while turning counter clockwise increases the gas and increases the flame.

You can adjust the flame temperature by rotating the AIR KNOB on the front of the unit. Rotating the knob reduces the air supply and lowers the flame temperature while turning counter clockwise increases the air supply and increases the flame temperature.

Connecting the foot switch

The footswitch is supplied with the StarFire ST

- Insert the jack of the footswitch in the socket on the back of the device (see page 7 for details)
- The foot switch is now connected and can be used to ignite the flame

Removing and attaching the burner head

- Pull the burner head up vertically to remove from base unit
- To reinstall, align the three contact pins with the proper holes and push gently straight down until it clicks in place

CAUTION

Do not lift or transport the StarFire ST by the burner head.

CAUTION

After using StarFire ST the burner head will be extremely hot, wait for this to cool down before removing the burner head.

Connecting StarFire ST to a power supply

- StarFire ST is supplied with a mains adapter power supply.
- Attach the mains adapter with the primary adapter suitable for your power line socket outlet.
- Plug in the mains adapter into a standard electrical source and connect it to the StarFire ST mains plug socket located in the back of the unit.

Operating StarFire ST

StarFire ST is on as long as there is a power supply.

Igniting the flame

StarFire ST can be ignited using the IGNITE button or the optional footswitch.

Using the ignite button

- To ignite the flame, briefly push the IGNITE button
- To extinguish the flame, briefly push the IGNITE button again

Using the footswitch

If the footswitch is plugged in (see 'Connecting the foot switch' on page 12) StarFire ST will automatically allow the use of the footswitch to ignite the flame.

- To ignite the flame, push the footswitch once
- The flame is extinguished upon release of the pedal

NOTE

While the footswitch is plugged in, the flame can be ignited or extinguished by either the IGNITE button or the footswitch.

Working with liquids

To minimize the risks of spilling liquids into or contaminating burner head, you may tilt the unit using the folding bracket located underneath the unit.

To tilt the device, pull the folding stand away from the base until it clicks into place. To un-tilt the device, push the folding stand together gently and rotate until flat against the base.

Maintenance, Care and Safety Precautions

Care and Maintenance

**WARNING**

Prior to cleaning make sure the unit is turned off and disconnected from the power source. Allow sufficient time for the burner head to cool. Disconnect the gas supply before proceeding.

StarFire ST can be cleaned with a soft cloth and most commercial disinfectants. A 70% ethanol solution can also be used.

The burner head can also be cleaned by following the removal procedure (see page 12) and cleaning with soap or commercial disinfectants.

Should the burner head fail to ignite you may try using a 70% ethanol solution applied with a nylon swab to clean the ignition electrode. Once the electrode is bent, however, it will no longer ignite and must be replaced.

**DANGER**

Before cleaning the equipment, remove the plug from the power supply. Never use aggressive solutions such as nitro thinners, etc.

Troubleshooting (FAQ)

In the event of an error the flame is automatically extinguished.

The following table is a list of the recommended action should StarFire ST extinguish the flame.

Problem	Remedy
Burner head does not spark or flame	<ul style="list-style-type: none"> • Check and make sure power is connected
Burner head sparks (audible click) but no flame	<ul style="list-style-type: none"> • Check to make sure gas is connected • If using gas cartridge ensure adapter has been turned to “on”. • Replace empty gas cartridge • Ensure the correct gas nozzle is in the unit • Verify burner head is not soiled or clogged • Inspect burner head and contact pins and clean when necessary. If electrode is bent or does not work after cleaning, replace burner head.
Flame blows out	<ul style="list-style-type: none"> • Adjust the gas/air mixture when using the controls on the front of the unit • Prevent excessive air flows or drafts • Check gas cartridge for fuel • Unit may have over heated due to continuous use. Allow the unit to cool for 30 minutes before attempting to restart.
Foot pedal does not work	<ul style="list-style-type: none"> • Disengage footswitch jack and then reinsert • Replace footswitch • Check to see if unit ignites by depressing the ignition button (see page 7)

If the problem still persists contact a service technician.

Technical Data

Equipment Type	Safety Bunsen Burner	
Dimensions (LxWxH)	178x122x39mm without burner head	
Weight	1.2Kg	
Power supply	Input voltage	100-240 VAC, 50/60Hz
	Output voltage	12VDC / 1.25A
Rechargeable Battery (optional)	Type	NiMH 3.6V
	Capacity	At least 2000mA
	Charging Time	12 h
	Operating Life	8 h
Operating Conditions	Operation	15-35°C, 15-80% RH (not condensing)
	Storage	-10-60°C, 5-80% RH (not condensing)
Nominal Thermal Rating	Propane/Butane gas	1.6kW
	Natural Gas	1.0kW
Nominal Gas Input Pressure	Propane/Butane gas	50mbar
	Natural gas	20mbar
Material	Top Casing	Stainless Steel
	Bottom Casing/ Battery Door	Zytel FR70M30
	Screen	Borosilicate Glass
	Feet	Rubber
	Knobs	Anodized Aluminium
Consumption	Propane/Butane gas	.12 Kg/h
	Natural Gas	.1 cbm/h
Rated Heat Load	Propane/Butane gas	1300W
	Natural Gas	1000W
Flame Temperature	Propane/Butane gas	1350°C
	Natural Gas	1300°C
CV360 Gas Cartridge Operation	40 min.	

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