MA160

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Vivaspin® Turbo 4 and 15

Allow fastest sample concentration
with highest recoveries

Benefits

- Complete Recoveries
- Fastest Concentrations
- Highest Chemical Resistance

Product Information

Vivaspin® Turbo 4 and 15 centrifugal concentrators offer the optimal solution to any concentration or buffer exchange application with their broad range of MWCOs.

Highest flow rates are achieved due to their twin vertical membranes which minimize protein polarization and subsequent fouling of the membrane. Additionally, their sleek internal profile ensures maximum process speeds right down to the last 100 µl.

Working Principle

Centrifugation provides the vector to clear solvent and micro molecules through an ultrafiltration membrane to separate macromolecular species and solvents primarily on the basis of size. It is particularly appropriate for the concentration of macromolecules and can also be used to purify molecular species or for solvent exchange. Ultrafiltration is a non denaturing method that is more efficient, flexible and gentle than alternative processes.

High Performance

In a single spin, 4 or 15 ml solutions can be concentrated up to 150x. Samples can be typically concentrated in 10–30 minutes with macromolecular recoveries in excess of 95%.

Optimized Design for Fastest Concentration Results

The Vivaspin® Turbo's optimized design, its sleek internal profile, ensures maximum process speeds right the way down to the last few micro liters. The UV joining technology allows for a smooth joint transition between membrane and plastic housing, allowing all of your valuable sample to be collected into the unique pipette friendly dead stop pocket.

The ultimate in centrifugal ultrafiltration technology:

– Sleek internal design

Ensures maximum process speed for the complete filtration

– Large twin vertical membranes

A fouling of the membrane is avoided due to minimized protein polarization.

– Unique angular dead stop pocket

The dead stop pocket in both Vivaspin® Turbo 4 & 15 is easy to access with standard 200 µl pipette tips due to its patent pending angular design. It eliminates the risk of the sample running to dryness while allowing highest recovery of the concentrate.

Optimized Choice of Materials for

High Chemical Compatibility

The combination of Polyethersulfone (PES), Polystyrene and Polypropylene (PP) allows sterilization and depyrogenization of the Vivaspin® Turbo 4 and 15 units.

Polyethersulfone membranes are preferred for their low fouling characteristics, exceptional flux and broad pH range compatibility.

Applications

Sample preparation

- Sample (protein, lipid, virus, nanoparticle, macromolecule) concentration
- Desalting | dialysis
- Buffer exchange

Your sample is often the result of several months of research. Your sample is valuable, and Vivaspin® Turbo 15 provides highest recovery.

Summary

For scientists and lab technicians who need to quickly and safely concentrate biological samples of 2 ml to 15 ml initial volume up to 150 fold, Sartorius offers the Vivaspin® Turbo 4 and 15 ultrafilters.

Unlike competitive ultrafiltration units, Vivaspin® Turbo 4 and 15 are equipped with an angular dead stop pocket, that enables reproducible and complete recoveries, while being the fastest in the market.

Technical specifications

		Vivaspin® Turbo 4	Vivaspin® Turbo 15
Materials	Body	Styrene butadiene copolymer	Styrene butadiene copolymer
	Filtrate vessel	Polypropylene	Polypropylene
	Concentrator cap	Polypropylene	Polypropylene
	Membrane	Polyethersulfone (PES)	Polyethersulfone (PES)
Dimensions	Total length (concentrator insert)	75.5 mm	77 mm
	Total length (in tube with cap)	122.5 mm	118 mm
	Diameter (concentrator insert)	14.6 mm	27 mm
	Active membrane area	3.2 cm ²	7.2 cm ²
	Hold up volume of membrane	<10 µl	<10 µl
	Dead stop volume in swing out	40 µl	100 µl
	Dead stop volume in fixed angle	30 µl	60 µl
Concentrator capacity	Swing bucket rotor	4 ml	15 ml
	Fixed angle rotor (25°)	4 ml	9 ml
Maximum speed 3 kDa – 50 kDa	Swing bucket rotor	4,000 × g	4,000 × g
	Fixed Angle rotor (25°)	7,500 × g	4,000 × g
Maximum speed 100 kDa	Swing bucket rotor	3,000 × g	3,000 × g
	Fixed Angle rotor (25°)	5,000 × g	3,000 × g
Sterilization	ETO or 70% EtOH		
Removal of endotoxins [Depyrogenization]	Flushing with 1N NaOH		

Performance Characteristics

Vivaspin® Turbo 4

Time to concentrate up to 30x [min.] at 20°C and solute recovery %				
Rotor	Swing bucket		Fixed angle (25°)	
Centrifugal speed	4,000 × g		7,500 × g	
Start volume	4 ml		4 ml	
	Min.	Rec.	Min.	Rec.
Cytochrome c* (12,400 MW)				
3 MWCO PES	60	98	80	96
Lysozyme* (14,300 MW)				
3 MWCO PES	65	95	70	93
Cytochrome c* (12,400 MW)				
5 MWCO PES	40	95	50	94
Lysozyme* (14,300 MW)				
5 MWCO PES	50	94	60	92
α-Chymotrypsin** (25,000 MW)				
10 MWCO PES	10	95	8	95
BSA** (66,000 MW)				
10 MWCO PES	10	98	7	97
30 MWCO PES	8	96	6	97
IgG** (160,000 MW)				
30 MWCO PES	18	94	13	92
50 MWCO PES	16	93	12	90
100 MWCO PES****	17	94	13	92

Vivaspin® Turbo 15

Time to concentrate up to 20x [min.] at 20°C and solute recovery %				
Rotor	Swing bucket		Fixed angle (25°)	
Centrifugal speed	4,000 × g		4,000 × g	
Start volume	15 ml		9 ml	
	Min.	Rec.	Min.	Rec.
Cytochrome c* (12,400 MW)				
3 MWCO PES	61	98%	86	97%
Lysozyme* (14,300 MW)				
3 MWCO PES	56	98%	87	97%
Cytochrome c* (12,400 MW)				
5 MWCO PES	30	98%	50	98%
Lysozyme* (14,300 MW)				
5 MWCO PES	33	96%	50	96%
α-Chymotrypsin** (25,000 MW)				
10 MWCO PES	10	95%	10	95%
BSA** (66,000 MW)				
10 MWCO PES	10	99%	10	99%
30 MWCO PES	8	98%	10	98%
IgG** (160,000 MW)				
30 MWCO PES	23	95%	17	95%
50 MWCO PES	20	94%	15	94%
100 MWCO PES***	30	92%	16	92%

* 0.25 mg/ml

** 1 mg/ml

*** 3,000 × g centrifugal speed

**** 3,000 × g swing-out | 5,000 × g fixed angle

Ordering Information

Vivaspin® Turbo 4 and 15 centrifugal concentrators, disposable ultrafiltration units, for processing sample volumes of 2 – 4 ml and 15 ml, Polyethersulfone UF membrane.

Vivaspin® Turbo 4

Cut off	Quantity	Order no.
3,000 MWCO	25	VS04T91
3,000 MWCO	100	VS04T92
5,000 MWCO	25	VS04T11
5,000 MWCO	100	VS04T12
10,000 MWCO	25	VS04T01
10,000 MWCO	100	VS04T02
30,000 MWCO	25	VS04T21
30,000 MWCO	100	VS04T22
50,000 MWCO	25	VS04T31
50,000 MWCO	100	VS04T32
100,000 MWCO	25	VS04T41
100,000 MWCO	100	VS04T42

Vivaspin® Turbo 15

Cut off	Quantity	Order no.
3,000 MWCO	12	VS15T91
3,000 MWCO	48	VS15T92
5,000 MWCO	12	VS15T11
5,000 MWCO	48	VS15T12
10,000 MWCO	12	VS15T01
10,000 MWCO	48	VS15T02
30,000 MWCO	12	VS15T21
30,000 MWCO	48	VS15T22
50,000 MWCO	12	VS15T31
50,000 MWCO	48	VS15T32
100,000 MWCO	12	VS15T41
100,000 MWCO	48	VS15T42

Sartorius Stedim Biotech GmbH
August-Spindler-Strasse 11
37079 Goettingen, Germany

Phone +49.551.308.0
Fax +49.551.308.3289
www.sartorius.com

USA Toll-Free +1.800.368.7178
UK +44.1372.737159
France +33.442.845600
Italy +39.055.63.40.41
Spain +34.90.2110935
Russian Federation +7.812.327.5.327
Japan +81.3.4331.4300

Specifications subject to change
without notice. Printed and copyrighted
by Sartorius Stedim Biotech GmbH. | W
Publication No.: SLU2018-e140606
Order No.: 85032-541-62
Ver. 06 | 2014

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www.sartorius.it/promo-optifit-tips

Sales and Service Contacts

For further contacts, visit www.sartorius.com

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Phone +3623.457.227
Fax +3623.457.147

Ireland

Sartorius Ireland Ltd.
Unit 41, The Business Centre
Stadium Business Park
Ballycoolin Road
Dublin 11

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Phone Spain +34.902.123.367
Phone Portugal +351.800.855.800
Fax Spain +34.91.358.96.23
Fax Portugal +351.800.855.799

Switzerland

Sartorius Mechatronics Switzerland AG
Ringstrasse 24a
8317 Tagelswangen (ZH)

Phone +41.44.746.50.00
Fax +41.44.746.50.50

U.K.

Sartorius UK Ltd.
Longmead Business Centre
Blenheim Road, Epsom
Surrey KT19 9QQ

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Asia | Pacific

Australia

Sartorius Australia Pty. Ltd.
Unit 5, 7-11 Rodeo Drive
Dandenong South Vic 3175

Phone +61.3.8762.1800
Fax +61.3.8762.1828

China

Sartorius Scientific
Instruments (Beijing) Co., Ltd.
33 Yu An Road, Airport Industrial Park
Zone B,
Shunyi District, Beijing 101300,
P.R.China

Phone +86.10.8042.6300
Fax +86.10.8042.6486

Hong Kong

Sartorius Hong Kong Ltd.
Unit 1012, Lu Plaza
2 Wing Yip Street
Kwung Tong
Kowloon, Hong Kong

Phone +852.2774.2678
Fax +852.2766.3526

India

Sartorius Weighing India Pvt. Ltd.
#69/2-69/3, NH 48, Jakkasandra,
Nelamangala Tq
562 123 Bangalore, India

Phone +91.80.4350.5250
Fax +91.80.4350.5253

Japan

Sartorius Stedim Japan K.K.
4th Fl., Daiwa Shinagawa North Bldg.
8-11, Kita-Shinagawa 1-chome
Shinagawa-ku, Tokyo, 140-0001 Japan

Phone +81.3.4331.4300
Fax +81.3.4331.4301

Malaysia

Sartorius Malaysia Sdn. Bhd
Lot L3-E-3B, Enterprise 4
Technology Park Malaysia
Bukit Jalil
57000 Kuala Lumpur, Malaysia

Phone +60.3.8996.0622
Fax +60.3.8996.0755

Singapore

Sartorius Singapore Pte. Ltd
1 Science Park Road,
The Capricorn, #05-08A,
Singapore Science Park II
Singapore 117528

Phone +65.6872.3966
Fax +65.6778.2494

South Korea

Sartorius Korea Ltd.
8th Floor, Solid Space B/D,
PanGyoYeok-Ro 220, BunDang-Gu
SeongNam-Si, GyeongGi-Do, 463-400

Phone +82.31.622.5700
Fax +82.31.622.5799

Thailand

Sartorius (Thailand) Co. Ltd.
129 Rama 9 Road,
Huaykwang
Bangkok 10310

Phone +66.2643.8361-6
Fax +66.2643.8367

sartorius stedim
biotech

Bioprocessing Media and Buffers Grow with Us

turning science **into solutions**

Experience

“ Highly experienced media on the selection of standard media as well as development This service packaging, documentation

Industry Leader

50 years of culture media experience backed by 150 years of pharma industry innovations

Transparency

Sharing experiences to bring added value and fit-for-purpose products

► Leading position in single-use applications including fluid management, filtration and process analytical technology

► Pioneer in state-of-the-art micronization high-quality powder production process

► Long-term relationships with dual-sourced raw material suppliers guarantees supply and quality
► Excellent regulatory dossier support

► Risk minimization approach applied to powder | liquid solutions
► ISO certification with successful audit track record following Current Good Manufacturing Practices 21CFR820

Florenz Sartorius mechanical workshop is founded

Sartorius starts commercial-scale manufacturing of membrane filters

Sartorius is founded as the original manufacturer of EVA plastic bags for total parental nutrition

Stedim creates the market for bag systems used in biopharmaceutical manufacturing processes

1870

1922

1948

1970

1974

1979

1990

1994

Microbiological associates


Whittaker Bioproducts: BioWhittaker Media

Grand opening of the Walkersville facility in Maryland, USA

Initial manufacture of liquid media in Verviers, Belgium, EU

Lonza

specialists will provide guidance
formulations,
and optimization of custom media.
includes product specific
and full regulatory support. ”

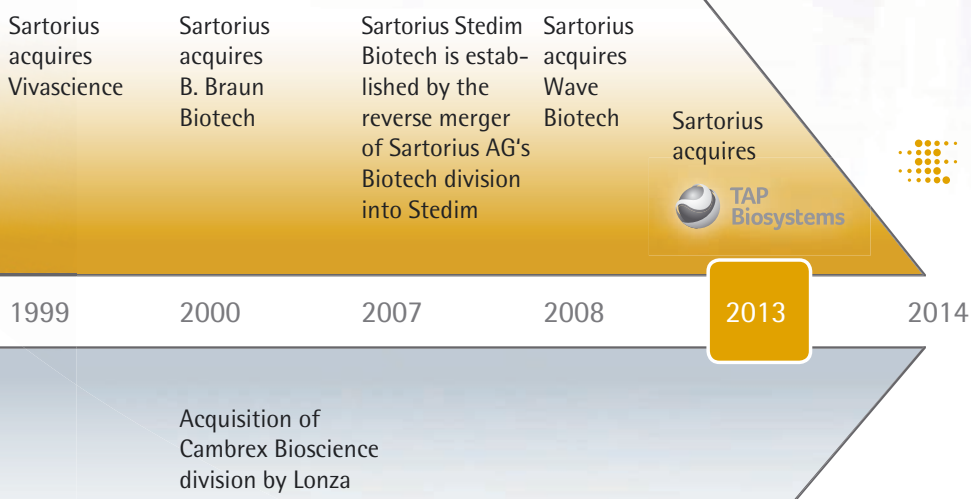Long-term Collaboration

Complete offering with
complementary product
lines to address all your
bioprocessing needs

Flexibility


Research

Development


Pre-Clinical

Phase I

Antibodies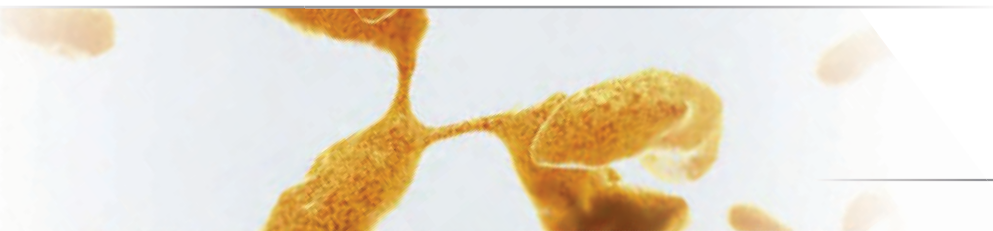


Proteins

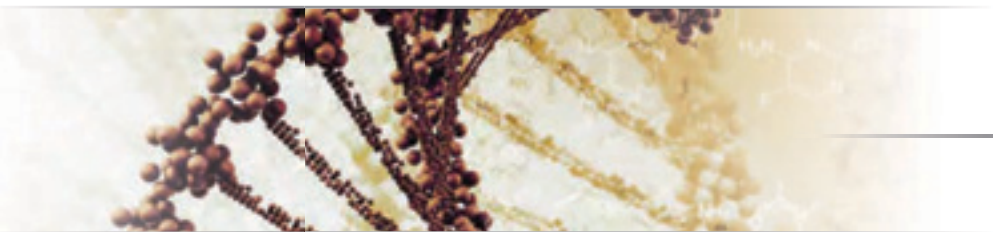Vaccines

Cell Therapy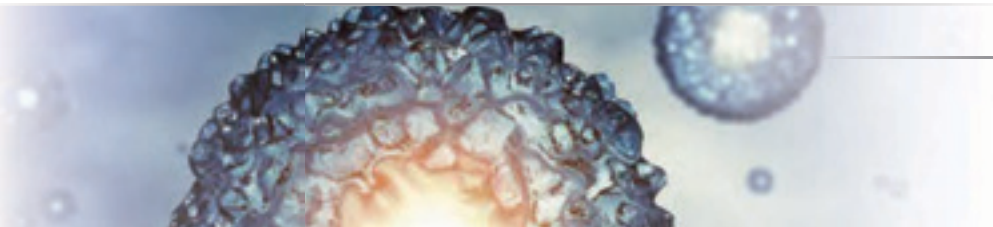

Hybridoma

CHO


Microorganisms

Insect Cells

Kidney Cell Line

Primary Cells

Stem Cells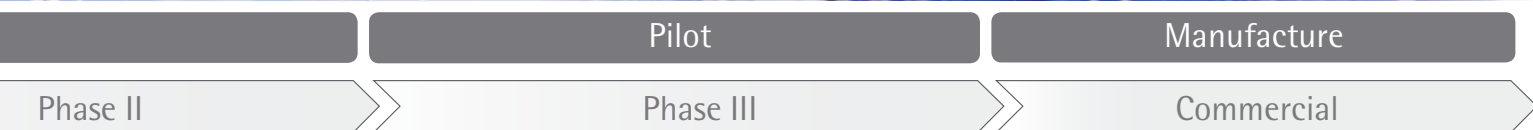


“ Working in collaboration, Lonza and Sartorius offer a wide array of liquid and powder media products along with the expertise to help scale up your project from clinical trials up to commercialization. Offered products and services include specially engineered media and buffers for dedicated cell lines and USP | DSP | F&F unit operations. ”

Industry leader in batch size scalability	▶ Liquid: 1 L to 10,000 L Powder: 7 kg to 7,000 kg
Dual global manufacturing sites	▶ Assurance of supply
Wide range of primary and secondary packaging	▶ For your specific needs
Off-the-shelf to fully customized solutions	▶ For maximized titers and full product life support
Formulation options include AO NAO*, Serum-free, protein-free with full traceability	▶ Solutions for any type of process

* AO | NAO : Animal Origin | Non-Animal Origin or Animal Derived Component Free

Powder

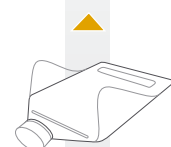5 to 20 kg

5 to 20 kg


100 g to 2 kg

10 g


Liquid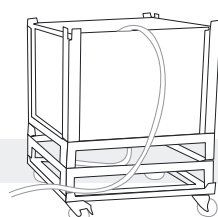

100 L to 1,000 L

50 L to 200 L

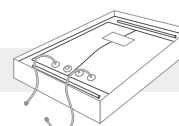1 L to 20 L

0.5 L to 2.5 L

Complete Solution

For cell line development and media selection ▶

Standards off-the-shelf media

High quality | high performance Serum free Formulation
for all cell lines in bottles and small bags

For media optimization ▶

Customized media optimization service and support

R&D Specialists can optimize or fully develop new custom media
engineered to your unique cell culture specifications

For process development ▶

Fast and clean custom production

An expedited production option for those that require high quality
cell culture media for process optimization or process development
as well as quick delivery of buffers

For production ▶

Custom media production under cGMP Protocols

Small to large batch production following Current Good manufac-
turing practices (cGMP 21CFR820) in ISO certified facilities
Manufacturing Practices compliance regulations with a standard
battery of product release tests on request

Processing Path Showing USP | DSP Elements

Media | Buffer Preparation

LevMixer

FlexAct®

Sterile Filtration

Storage

Biostat® B, RM, UniVessel

Upstream Process

Pre-Clinical

Media

Pre-Viral Downstream Process

Protein A Chromatography

Low-pH Virus Inactivation

CEX Chromatography

Storage

Wash | Equilibration | Elution buffers

pH Adjustments buffers

Wash | Equilibration | Elution buffers

Post-Viral Downstream Process

FlexAct®

UF | DF

Storage

Wash | Flush | Sanitize | Diafiltration buffers

A Media Solution to Fit Your Cell Line, Schedule and Budget

Features

Media and buffers in Sartorius Stedim bags

Full media qualification package and support. All liquid media filtered through Sartorius filters

Prew weighed powder buffer and media in dispensing bags

Integrated media feed strategies and PAT automation

Products fully integrated into the Sartorius upstream and downstream solutions

Benefits

- ▶ Same material as for single-use bioreactor bags to minimize the risk of leachables
- ▶ Optimized media filtration and guaranteed successful scale-up thanks to complete support provided
- ▶ Simple, ready-to-use media and buffers fully compatible with mixing technology
- ▶ Fully monitored and controlled cell culture conditions
- ▶ One-stop supplier gives you full support so you can focus on your core tasks

➤ Cell Growth Media ➤ Buffers

Biostat® STR

FlexAct®

Depth Filtration

Polishing

Storage

Feeds

Wash buffers

Wash | Equilibration | Elution buffers

Polishing

Storage

FlexAct®

Virus Filtration

Storage

Flush | Equilibration buffers

Buffer Flush buffers

Sterile Filtration

Filling

Freezing

Dilution | Stabilization buffers

Sartorius Stedim Biotech and Lonza Global Facilities

For more information about our collaboration and for a complete list of our offices worldwide, please visit www.sartorius-stedim.com/media-buffer

Sartorius Stedim Biotech Facilities

Goettingen

August-Spindler-Strasse 11, 37079 Goettingen, Germany
Phone +49.551.308.0 | Fax +49.551.308.3289

Guxhagen

Robert-Bosch-Strasse 5-7, 34302 Guxhagen, Germany
Phone +49.5665.407.0 | Fax +49.5665.407.2200

Aubagne

ZI des Paluds, Avenue de Jouques – CS 91051
13781 Aubagne Cedex, France
Phone +33.442.845600 | Fax +33.442.845619

New York

5 Orville Drive, Bohemia, NY 11716, USA
Toll-Free +1.800.368.7178 | Fax +1.631.254.4253

Lonza Facilities

Lonza

Biotherapeutic Cell
Culture Media

Verviers

Parc Industriel de Petit-Rechain, BE-4800 Verviers, Belgium
Phone +32.8732.1611 | Fax +32.8735.1967

Walkersville

8830 Biggs Ford Road, Walkersville, MD 21793, USA
Phone +1.301.898.7025 | Fax +1.301.845.4024

◀ www.sartorius-stedim.com/media-buffer

Packaged WFI (Water for Injection) Products

Our Water For Injection (WFI) products are manufactured and tested according to strict industry standards. It meets or exceeds USP and EP requirements.

2014 Promotion 20 L and 200 L Bags

- Produced according to Current Good Manufacturing Practices conditions
- US and European production plants act as back up with similar technologies
- Can be customised (larger volumes, 3D bags)
- Consistent Production Methods and Documentation for all sizes
- Complementary
 - For mixing solution
 - Filter products (washing)

Definition

Our off-the-shelf WFI products are available in 500 ml or 1 L bottles, 20 L and 200 L bags. We can also provide made to-order WFI in custom packaging options to fit your application or direct ready-to-use solutions (Cell Culture Media, Buffers, Organ Transplant Solutions).

Delivery time: 24 hours to 1 week from reception of Purchase Order (Add an extra week for Asia).

Facilities follow cGMP 21CFR820 and are certified ISO9001 and ISO13485.

Product Description

Water For Injection produced in bulk; meeting USP and EP Pharmacopeia criteria; filtered using Sartopore II filters into primary packaging (bottles and | or bags)

– Available in 500 mL or 1 L bottles, in 20 L or 200 L Bags off the shelf

Bulk packaged for commercial use elsewhere, not for parenteral administration. Our WFI Quality Water is aseptically processed from a validated Water For Injection System that meets current USP and EP specifications in the loop. Our WFI Quality Water is not to be confused with Sterile Water for Injection which is intended for use in extemporaneous prescription compounding, and as a diluent for injectable parenteral products. In addition, Sterile Water for Injection must be packaged in single dose containers not larger than 1 liter in size.

This product is manufactured aseptically under cGMP 21CFR820 conditions by a validated sterile filtration method and tested according to current U.S.P. and E.P.

This product has been produced for laboratory | manufacturing use. Use for in vitro diagnostic procedures has not been established. It is the end user's responsibility to ensure that the final products meets the requirements of the application.

Application

Subject to the proviso, customer checks quality according to CoA is suitable for his needs

UpStream: Media Preparation and Hydration


Downstream: Buffer Preparation

UpStream & Downstream: Cleaning, Washing, Rinsing (filters, equipment)

Features	Benefits
Very Low Endotoxin level and Sterile	Minimization of Contamination
High Quality made under cGMP 21CFR820 conditions	Regulatory Friendly and Risk Minimization
No need for internal installation of WFI loop	Significant Savings

Certificate of Analysis (CoA)

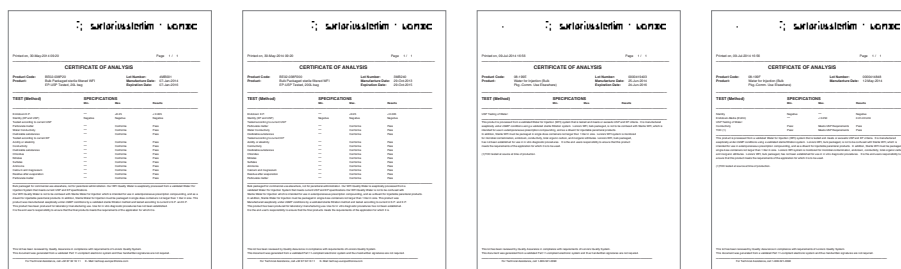CoA included with all product sold. CoA is region and packaging size dependent.

Ordering Information

Part No.	Description
BE17-724F	500 mL Bottle, Water for Cell Culture-WFI Quality
BE17-724Q	1 L Bottle, Water for Cell Culture-WFI Quality
BE02-038P20	20 L Bag, Water for Cell Culture-WFI Quality
BE02-038P200	200 L Bag, Water for Cell Culture-WFI Quality

Sartorius Stedim Biotech GmbH
August-Spindler-Strasse 11
37079 Goettingen, Germany

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* Tablet | smartphone non inclusi

Sales and Service Contacts

For further contacts, visit www.sartorius-stedim.com

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Phone +41.52.354.36.36
Fax +41.52.354.36.46

U.K.

Sartorius Stedim UK Ltd.
Longmead Business Centre
Blenheim Road, Epsom
Surrey KT19 9 QQ

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Phone +61.3.8762.1800
Fax +61.3.8762.1828

China

Sartorius Stedim Biotech (Beijing) Co. Ltd.
No. 33 Yu'an Road
Airport Industrial Park Zone B
Shunyi District, Beijing 101300

Phone +86.10.80426516
Fax +86.10.80426580

Sartorius Stedim Biotech (Beijing) Co. Ltd.
Shanghai Branch Office
3rd Floor, North Wing, Tower 1
No. 4560 Jin Ke Road
Pudong District, Shanghai 201210

Phone +86.21.68782300
Fax +86.21.68782332 | 68782882

Sartorius Stedim Biotech (Beijing) Co. Ltd.
Guangzhou Representative Office
Unit K, Building 23
Huihua Commerce & Trade Building
No. 80 Xianlie Middle Road
Guangzhou 510070

Phone +86.20.37618687 | 37618651
Fax +86.20.37619051

India

Sartorius Stedim India Pvt. Ltd.
#69/2-69/3, NH 48, Jakkasandra
Nelamangala Tq
562 123 Bangalore, India

Phone +91.80.4350.5250
Fax +91.80.4350.5253

Japan

Sartorius Stedim Japan K.K.
4th Fl., Daiwa Shinagawa North Bldg.
8-11, Kita-Shinagawa 1-chome
Shinagawa-ku, Tokyo, 140-0001 Japan

Phone +81.3.4331.4300
Fax +81.3.4331.4301

Malaysia

Sartorius Stedim Malaysia Sdn. Bhd.
Lot L3-E-3B, Enterprise 4
Technology Park Malaysia
Bukit Jalil
57000 Kuala Lumpur, Malaysia

Phone +60.3.8996.0622
Fax +60.3.8996.0755

Singapore

Sartorius Stedim Singapore Pte. Ltd.
1 Science Park Road,
The Capricorn, #05-08A,
Singapore Science Park II
Singapore 117528

Phone +65.6872.3966
Fax +65.6778.2494

South Korea

Sartorius Korea Biotech Co., Ltd.
8th Floor, Solid Space B/D,
PanGyoYeok-Ro 220, BunDang-Gu
SeongNam-Si, GyeongGi-Do, 463-400

Phone +82.31.622.5700
Fax +82.31.622.5799