Agitador Vórtex 681/2 | Vortex mixer 681/2

Marca|Trademark: Nahita

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Nahita vortex mixer 681/2 allows a quick and effective mixture of substances contained in a wide range of recipients, thanks to its orbital vibrating movement. A special anti-sliding four feet system and its flat construction ensure a high lab bench stability while using the instrument, and avoid transmission of vibrations to other near equipments.

Speed can be adjusted to up to 1800 rpm for a best fit to end-user requirements. Nahita vortex mixer includes also a double operational working mode, continuous and discontinuous-touch mode (mixing activity starts by a light pressure of the test tube on the attachment), and is equipped with different accessories easily interchangeable that give the equipment a high flexibility of use:

- Attachment for test tubes up to 30 mm diameter. A trough prevents slipping caused by vibrations. For using in touch mode.
- Foam attachment for 6 Eppendorf-type tubes (1.5 mL). For using in continuous mode.
- Attachment for microtiter plates. For using in continuous mode.
- Foam attachment for Petri dishes, culture flasks and Erlenmeyers. For using at a low shaking rate in continuous or discontinuous mode.
- Foam attachment for 6 test tubes of 20 mm diameter. For using in continuous mode.

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This manual should be available for all users of these equipments. To get the best results and a higher duration of this equipment it is advisable to read carefully this manual and follow the processes of use.

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INDICE DE IDIOMAS

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5. MAINTENANCE AND CLEANING

To get the best results and a higher duration of this equipment it is essential to follow the processes of use.

Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- ♦ Please follow the processes of use of this manual.
- ♦ This manual should be available for all users of this equipment.
- ♦ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ♦ The vortex mixer is supplied with a Schuko standard wire. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ♦ Never try to repair the vortex mixer by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the vortex mixer (burns, hurts...) or damages in nearby equipments.
- ♦ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

Cleaning

- ♦ Never use scratching scourers or substances for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the vortex mixer and produce an early ageing of the equipment.
- ♦ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

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Security

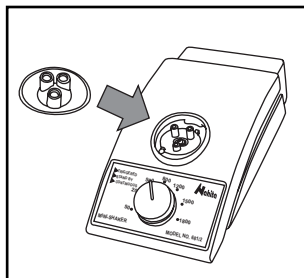
- ♦ The vortex mixer must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ♦ Please follow the relevant safety instructions and guidelines and occupational health and safety regulations for use in the laboratory
- ♦ When working with the shaking unit, the user must select and wear his personal protective equipment according to the hazard category of the mixing medium. Defective or inappropriate protective equipment can expose the user to the risk of spurting liquids, projectile parts or being pulled in at the shaking table or support.
- ♦ You should put the vortex mixer in a horizontal plane stable table, having a safety area of at least 30 cm per side.
- ♦ The feet of the appliance must be clean and undamaged
- ♦ Before use, please ensure that the knob for setting the speed is at the left stop as the appliance starts up at the last speed which was set.
- ♦ Ensure that parts of the body (hair or items of clothing) cannot be trapped by the moving parts.
- ♦ All accessories and vessels in place for the shaking process must be firmly secured. Shaking vessels which are not properly secured could get damaged or be projected out, thus causing injury. It is essential to regularly check that the vessels to be shaken and the attachments are firmly secured, especially before using the appliance again.
- ♦ If you notice that the device is not running smoothly, the speed must always be reduced until no more unevenness occurs in the operation. Please make certain that individual agitation tubes are placed in the middle of the agitation table and that if several tubes are placed on the agitation table together, they are evenly distributed and well secured.
- ♦ During its functioning dangerous materials such as flammable or pathological substances must be out of the safety area.
- ♦ The appliance starts up again automatically following a cut in the power supply during a work process.
- ♦ When you are not using the vortex mixer for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ♦ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ♦ Never try to repair the vortex mixer by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ♦ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the vortex mixer (see Maintenance).
- ♦ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

ATTENTION! Before use, please ensure that the knob for setting the speed is at the left stop as the appliance starts up at the last speed which was set.

Placing attachments

- ♦ As a standard, the vortex mixer includes an adapter for 6 tubes of 20 mm diameter and several attachments: for 6 tubes Eppendorf-like, for tubes up to 30 mm diameter, for microtiter plates and an attachment for flasks.
- ♦ Insert the desired attachment, fitting the three projections that it has with the holes of the equipment body. Make sure the attachment is well fitted.
- ♦ Please insert the wire that feeds the A/C to the base of current 220V $\pm 10\%$ 50-60Hz provided with earth wire and, on the other extreme, to the mixer vortex. Switch the mixer vortex. If you want to work in continuous operation mode, set the switch in position (I) and if you want to work in touch mode, set the switch in position (II).


Continuous operation mode (I)

a) Working with attachment for tubes up to 30 mm:

- ♦ To set the vortex mixer to continuous mode, set the switch in position (I) and hold with the hand the tube on the platform. Adjust the desired speed.
- b) Working with attachment for tubes of 20 mm:*
- ♦ Put the tubes into the attachment, switch the mixer vortex and adjust the speed.

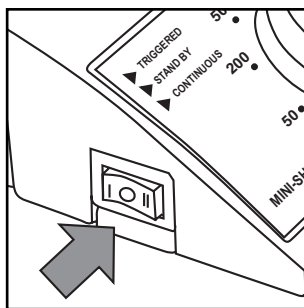Note: You should secure the accessories and vessels firmly in place, otherwise shaking vessels could be damaged.

Touch mode

- ♦ To set the vortex mixer to touch mode, set the switch in position (II)
- ♦ Adjust the desired speed with the speed knob (4) and press a tube gently with the hand on the platform.
- ♦ This way of working is suitable for shaking in press mode without changing the shaking speed, which enables making more than one in few time.
- ♦ Once the shaking is finished turn off the mixer vortex, switch in position (0) and unplug it.

Note: Reduce the speed if:

- The medium splashes out of the vessel because the speed is too high.
- The appliance is not running smoothly.
- The appliance begins to move around because of dynamic forces.

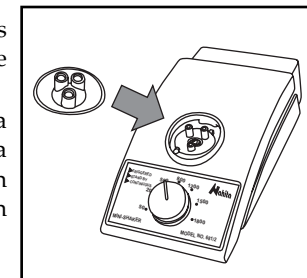
- ♦ If you are not using the vortex mixer for a long period of time please make sure it is disconnected from the net and protected from dust (this way you will avoid accidents and will extend its working-life).

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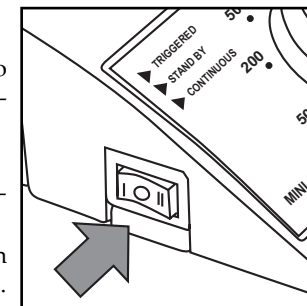
4. INSTALLATION / SETTING UP**Preliminary inspection**

- ♦ Unwrap the vortex mixer, take off the involving plastic and take off the polispán protection in which it comes fitted. Take off all the protective items and, without connecting the vortex mixer to the net, make sure that it does not present any damage because of the shipment. In case the equipment presents any damage tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.
- ♦ Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.
- ♦ Please check that all the accessories are enclosed with the equipment:
 - Vortex mixer
 - Adapter for 6 tubes of 20 mm diameter
 - Attachment for 6 tubes Eppendorf-like.
 - Attachment for tubes up to 30 mm diameter
 - Attachment for microtiter plates
 - Attachment for flasks
 - Shuko wire
 - User manual
 - Warranty

We will not accept any equipment in return period unless it comes in its original wrapping.

Installation

- ♦ Before using this instrument, it is convenient for you to familiarize with its components and basic essentials.

PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT.

- ♦ Please put the vortex mixer on top of a horizontal, plane and stable table making a free space of at least 30 cm per side. Do not put the vortex mixer near any warm supply (burners, blowlamps...), nor expose it to direct sunlight.
- ♦ Avoid inflammable or toxic substances in the working area.
- ♦ The vortex mixer is supplied with a Schuko standard wire
- ♦ Please insert the wire that feeds the AC electric current in the base of current 220V 50Hz $\pm 10\%$ provided with earth wire and to the other end to the vortex mixer connector.

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to people suffered by the improper use of the electric connection. The tension should be 220V 50Hz $\pm 10\%$.

- ♦ With Nahita vortex mixer you can chose continuous and discontinuous shaking mode as you need.


INDEX OF CONTENTS

1. USES OF THE INSTRUMENT	10
2. DESCRIPTION	10
3. TECHNICAL SPECIFICATIONS	10
4. INSTALLATION / SETTING UP	11
5. MAINTENANCE AND CLEANING	14
ANNEX I: CE CERTIFICATE	15

1. USES OF THE INSTRUMENT

This little sitrrer gets a fast mixture of the substances contained in a test tube, thanks to its vibratory movement. It is designed for use in laboratories, chemical industry, schools, pharmacies.

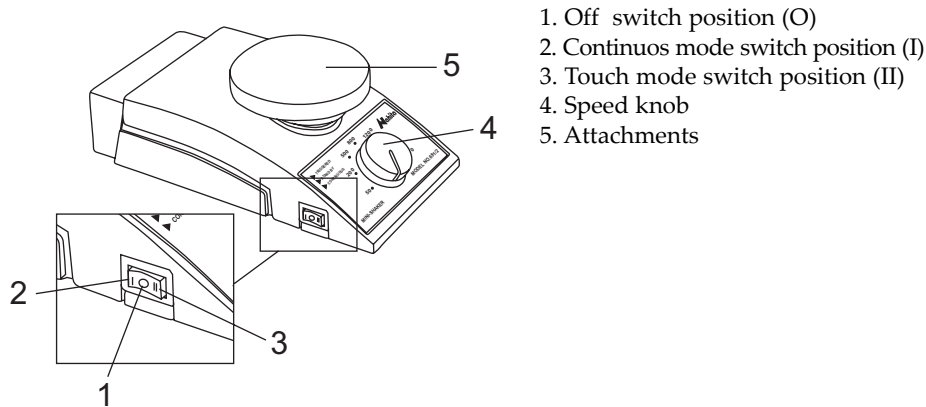2. DESCRIPTION

3. TECHNICAL SPECIFICATIONS

Code	50681002
Continuous mode	Yes
Touch mode	Yes
Speed range	50-1800 r.p.m

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www.twistermedical.com
info@twistermedical.com
902 876 983

Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita vortex mixer. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user. The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of *Maintenance*, where we expose the security notes.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Please bear in mind the following:

- ♦ This manual is inseparable from the nahita vortex mixer, so it should be available for all the users of this equipment.
- ♦ You should carefully handle the vortex mixer avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ♦ Never dismantle the different pieces of the vortex mixer to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the vortex mixer.
- ♦ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ♦ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to asistencia@auxilab.es or by telephone: +34 807 117 040 (0,30 Euros/min).
- ♦ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ♦ Overhaul is not covered by the vortex mixer warranty.
- ♦ Operations made by non-qualified staff will automatically produce a loss of the vortex mixer warranty.
- ♦ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ♦ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ♦ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.
- ♦ Manufacturer reserves the right to modify or improve the manual or equipment.

ATTENTION!

IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

INDEX OF LANGUAGES

Spanish	2-8
English	9-15