**ROTAVAL
80040**

EVAPORADOR ROTATORIO AL VACÍO

VACUUM ROTARY EVAPORATOR

ROTAVAL

MANUAL

Indice/Contents

INFORMACIÓN GENERAL	4
GENERAL INFORMATION	4
LISTA DE EMBALAJE	4
PACKING LIST	4
ESPECIFICACIÓN TÉCNICA.....	6
TECHNICAL FEATURES	6
INSTALACIÓN	6
INSTALLATION.....	6
OPERACIÓN	8
OPERATION.....	8
MANTENIMIENTO	11
MAINTENANCE	11
GARANTÍA.....	11
GUARANTEE	11
DECLARACIÓN DE CONFORMIDAD “CE”	12
“EC” CONFORMITY DECLARATION	12

INFORMACIÓN GENERAL

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or the ICC technical service.*
- 5) **IMPORTANT! ICC WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*

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PACKING LIST

The standard equipment consists of the following components:

Rotation and vacuum unit	
Heating bath	
Cooling bath with pump	
Temperature probe condensator	
Instruction manual	80214

EQUIPMENT PARTS:

35	2-litre evaporation bottle
36	1-litre collecting bottle
37	Screwdriver
38	Joint gasket
39	Fixing grips
40	Fuses set
41	Power supply cord
42	Heating bath cord
43	1-litre collecting bottle
44	1-litre evaporation bottle
45	3-litre collecting bottle
46	Valve
47	Adapter to small evaporation bottles
48	Condensator

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TECHNICAL FEATURES

<i>Voltage supply</i>	230V 50 Hz
<i>Volume of evaporator</i>	50-3000 ml
<i>Rotor rack up and down range</i>	0-150 mm
<i>Rotary speed</i>	20-200 rpm
<i>Rotary power</i>	40 W
<i>Bath heating tank measurement:</i>	
<i>Height</i>	14 cm
<i>Diameter</i>	28 cm
<i>Heating bath power</i>	1400 W
<i>Heating bath temperature range</i>	+5 ~ 90°C
<i>Overall dimensions:</i>	
<i>Height</i>	79 cm
<i>Width</i>	58 cm
<i>Depth</i>	45 cm
<i>Weight:</i>	37.2 kg

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INSTALLATION

Place the equipment on a flat, horizontal and level surface.

Assemble the supplied tubes and flasks.

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Be sure that the voltage supply is the same as the one indicated on the characteristic plate of the apparatus.

Do not use the apparatus if it is NOT earthed.

Unplugg the equipment from the power supply before replacing a fuse.

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CONNEXIONS AND PIPING:

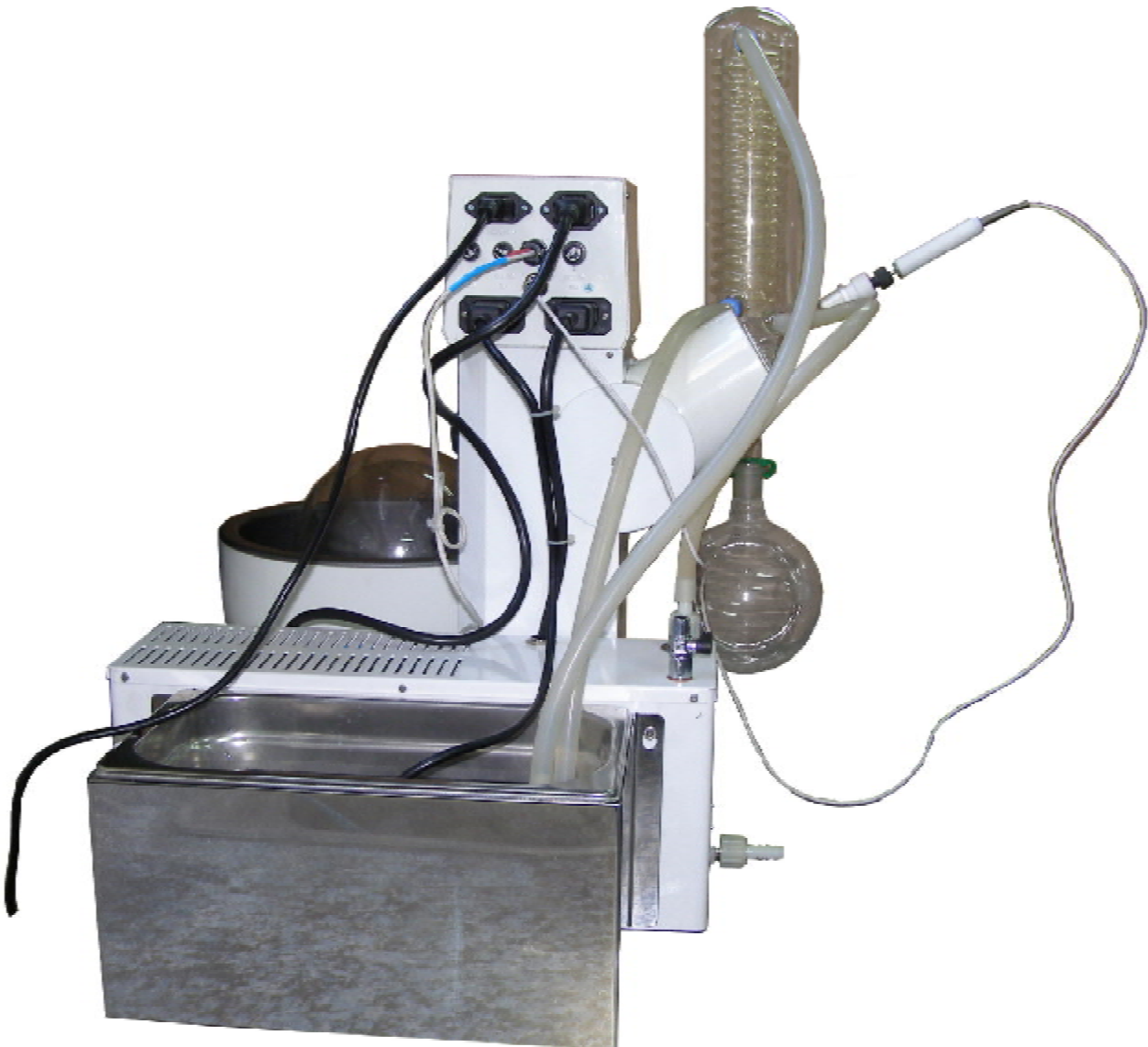
Connect the equipment to the power supply. Connect the heating bath and the water circulation pump, as it is showed on the picture, taking as a guidance the descriptions in the section OPERATION.

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IT IS ADVISABLE TO KEEP THE UNION BETWEEN THE GLASSES AND THE ROTOR CLEAN AND GREASED WITH FAT OR OIL COMPATIBLE WITH THE FOOD USE

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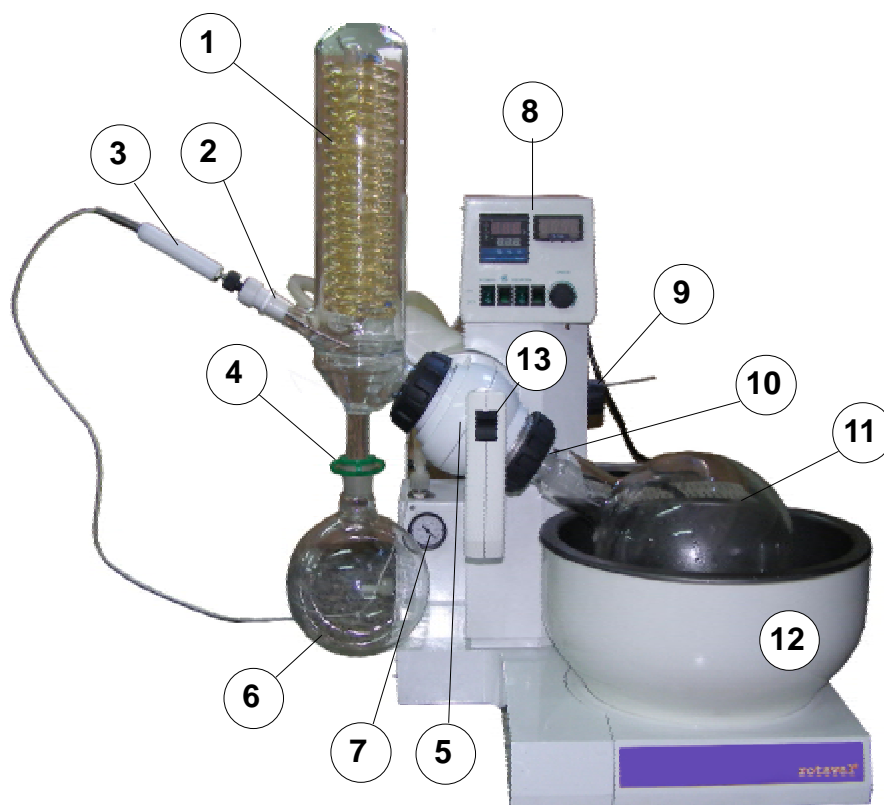
OPERATION

CAUTION BEFORE USING:

- The user must read this instruction manual
- When you need to work with an expensive product, it is advisable to try before with a small sample.
- The water temperature of the bath is controlled automatically and showed on the digital display.
- The evaporator must not be used in flammable ambients.

EQUIPMENT PARTS:

- 1 Condensator
- 2 Probe inlet
- 3 Condenser temperature probe
- 4 Fixing grip
- 5 Rotor
- 6 Collecting bottle
- 7 Vacuum meter
- 8 Control panel
- 9 Inclined angle adjustor
- 10 Locking knob
- 11 Evaporating flask
- 12 Water heater
- 13 Elevator push button

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
REAR PART

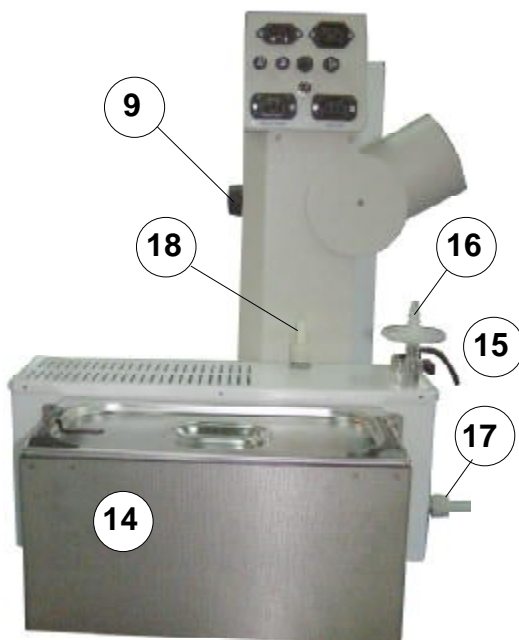
- 14 Cold bath
- 15 Vacuum regulation valve
- 16 Air inlet filter
- 17 Vacuum pump air expulsion
- 18 Vacuum pump aspiration

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CONTROL PANEL:

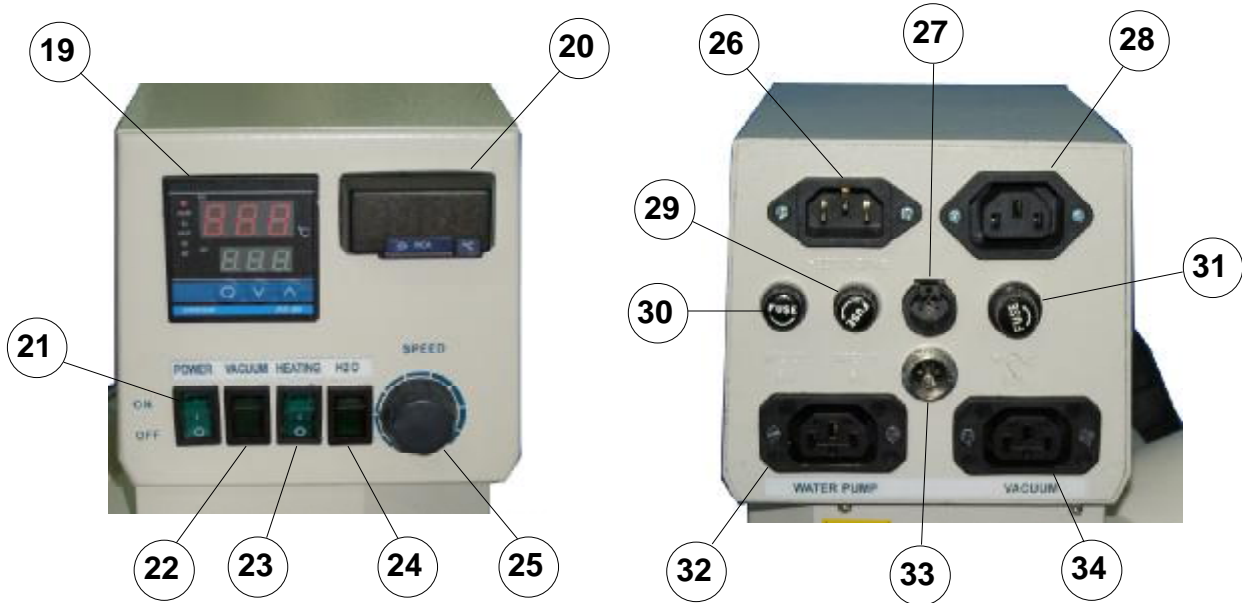- 19 Bath temperature control
- 20 Condensator temperature indicator.
- 21 Main switch
- 22 Vacuum pump switch
- 23 Heating bath swith
- 24 Condensator water circulation swith
- 25 Rotor speed regulatin knob
- 26 Mains connector
- 27 Heating bath probe connector
- 28 Heating bath connector
- 29 Fuse
- 30 Fuse
- 31 Fuse
- 32 Water pump connector
- 33 Condensator pump connector
- 34 Vacuum pump connector

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STARTING UP:

1. Be sure that the pipes are connected, temperature probe of the condenser (3) placed and the inlet air filter (16) inserted.
2. Fill water in the heating bath (12).
3. Turn on the main switch (21)
4. Set the working temperature on the temperature controler (19).
5. Turn on the switch (23) to start the heating
6. Fill in the cooling bath (14) with water and ice or antifreezing liquid.
7. Fill the liquid to be evaporated in the bottle (11)
8. Fix the bottle (11) with the nut (10)
9. Place the bottle (6) and loock it with the grip (4)
10. Adjust the position of the bottle (11) with the knob (9) in the way that do easier its contacts with the with the water of the heating bath (12).
11. Make lower the evaporation bottle (11) with the push button (13) untill the evaporation bottle be is slightly in the water of the heating bath (12).
12. Turn on the vacuum switch (22) and the water circulation (24).
13. Adjust the vacuum level with the manual valve (15) and rotation speed with the knob (25)

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MAINTENANCE

Before removing the cover disconnect the apparatus from the mains.

The control elements and heating elements must only be manipulated by qualified personnel.

CLEANING:

Keep clean and greased the close joints of the conical conical ground glass point with the condenser and the evaporation bottle. Keep the glasses cleaned.

GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use of causes beyond the control of ICC..

Any manipulation of the apparatus by unauthorized personnel by ICC. cancels the guarantee automatically.

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