**CENTRÍFUGA MEDIBAS
MEDIBAS CENTRIFUGE
CENTRIFUGEUSE MEDIBAS**

MODELO 2740 / MODEL 2740 / MODÈLE 2740

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**INSTRUCTIONS ON ENVIRONMENT PROTECTION**

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

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Thank you for choosing this equipment. We sincerely wish that you enjoy your centrifuge. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Please bear in mind the following:

- ◆ This manual is inseparable from the Nahita Medibas 2740 centrifuge, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the centrifuge avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the centrifuge to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the centrifuge.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to asistencia@auxilab.es
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the centrifuge warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the centrifuge warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

7. TROUBLESHOOTING

TROUBLE	CAUSE	SOLUTION
Displays do not turn on or turn on not correctly	<ol style="list-style-type: none"> 1. Fuse blown 2. Problem with the connection cable of the PCB 3. Problems with the chip of the equipment 	<ol style="list-style-type: none"> 1. Change the fuse 2. Contact with Nahita Technical Service 3. Contact with Nahita Technical Service
Displays turn on but equipment cannot operate	<ol style="list-style-type: none"> 1. Lid is not well closed 2. After the last operation, the lid has not been open. 3. After pressing Run, the button does not make good contact. 	<ol style="list-style-type: none"> 1. Close properly the lid 2. Open the lid, close it properly and try again. 3. Contact with Nahita Technical Service
The centrifuge makes too much noise	<ol style="list-style-type: none"> 1. The equipment is not well leveled 2. The surface in which the equipment is placed is not stable. 3. Mechanical components loose 4. Motor damaged 5. Any of the adapters does not swing out properly 6. Tubes have not been placed completely balanced 7. Dirt present on the rotor or inside the adapters 	<ol style="list-style-type: none"> 1. Put the equipment on top of a flat surface and level it correctly. 2. Put the equipment on top of a flat and stable surface 3. Contact with Nahita Technical Service 4. Contact with Nahita Technical Service 5. Check that all the adapters are correctly placed and move freely 6. Place again tubes in the adapter making sure they are perfectly balanced. 7. Clean properly rotor and adapters
Unstable rotating speed	<ol style="list-style-type: none"> 1. Failure in the control circuit 2. Brushes are worn out or copper bars are dirty 	<ol style="list-style-type: none"> 1. Contact with Nahita Technical Service 2. Contact with Nahita Technical Service

- ◆ Tube adapters are made of Nylon PA66 and can be sterilized in autoclave.
- ◆ Remember that the adapter buckets supplied together with the centrifuge wear out with the use; in case this happens, it is recommended to replace all adapters at the same time to avoid rotor unbalance.
- ◆ Never force the stop of the centrifuge, it is a risk of accident and debilitation and it may produce as well a mixing of the separated parts.
- ◆ Change of fuse: in case it is necessary to replace the fuse, it is situated near the socket.
- ◆ The change of brushes must be done by the Nahita Technical Service when the length is less than 6 mm.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ◆ The centrifuge is supplied with a Schuko standard wire. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ◆ Never try to repair the centrifuge by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the centrifuge (burns, hurts...) or damages in nearby equipments.
- ◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

Cleaning

- ◆ When working at high speed there is a risk of rupture of the tubes. If this happens, the stainless steel centrifuge chamber, the adapters and the buckets must be cleaned and disinfected with a disinfectant agent and the broken parts must be replaced; otherwise different components of the centrifuge will suffer important damages and due to the lack of cleaning mistakes in subsequent analyses will take place.
- ◆ For the cleaning of the different parts of the equipment, we recommend the use of neutral products with a pH value not higher than 6. These parts are also autoclavable.
- ◆ Pay special attention on aluminium alloy components; periodic treatment with antioxidant agents will prolong its life-span.
- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the centrifuge and produce an early ageing of the equipment.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

INDEX OF LANGUAGES

Spanish	2-12
English	13-23
French	24-34

INDEX OF CONTENTS

1. USES OF THE EQUIPMENT.....	14
2. DESCRIPTION	15
3. TECHNICAL SPECIFICATIONS	16
4. THEORETICAL FUNDAMENTALS.....	16
5. INSTALLATION / SETTING UP	17
6. MAINTENANCE AND CLEANING	20
7. TROUBLESHOOTING	22
8. ADAPTABLE ACCESSORIES RECOMMENDED	23


1. USES OF THE INSTRUMENTSCRIPCIÓN

The Nahita Medibas 2740 centrifuge is a bench-top, silent and sturdy model without vibrations. It is equipped with a unique swing-out rotor and three types of swinging adapters with capacities for 4 tubes of 100 ml (100x40 ml), 8 falcon standard-like tubes of 50 ml and 16 falcon standard-like tubes of 15 ml. The maximum speed the centrifuge can reach is 4000 r.p.m.

Because of its simple conception and the quality of the materials this centrifuge is specially indicated for any kind of analysis in Biology, Chemistry, Biochemistry, Veterinary... It can remain operating for long periods of time or keep a high daily rate of cycles without any difficulty.


2. DESCRIPTION

- 1.1 Cover
- 1.2 Rotor
- 1.3 Control panel
- 1.4 Electric current
- 1.5 Fuse
- 1.6 General switch ON/OFF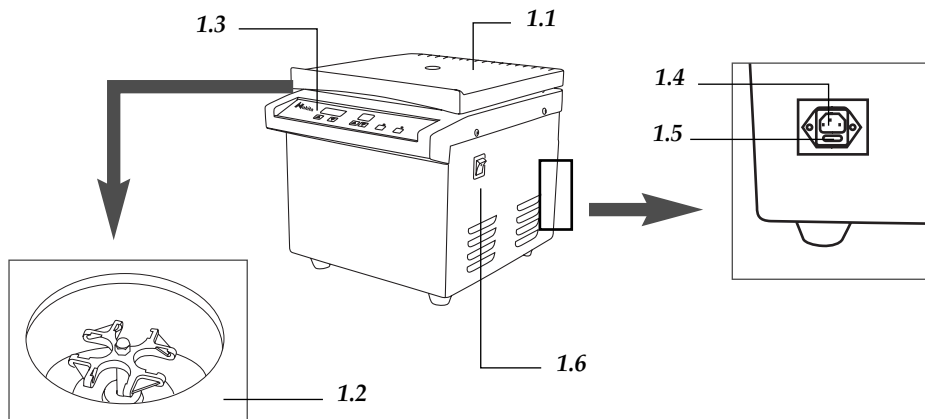

Front panel

- 2.1 Display of speed ("Speed")
- 2.2 Selector increase / decrease speed ▼▲
- 2.3 Display of time ("Time")
- 2.4 Selector increase / decrease time ▼▲
- 2.5 Run: start button
- 2.6 Stop: stop button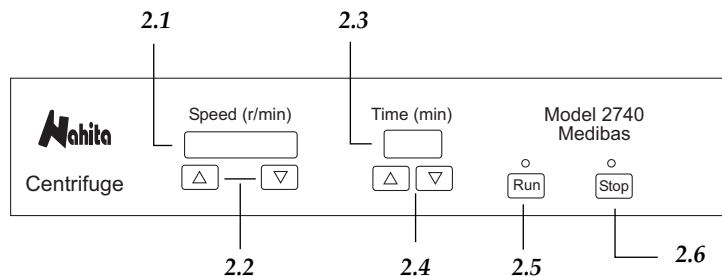

Security

- ◆ The centrifuge must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ Since it is very heavy, take the necessary precautions to assure the equipment is properly handled.
- ◆ During operation, dangerous material such as flammable or infectious liquids must be out of the working area.
- ◆ Nahita centrifuges are not suitable for treatment of flammable liquids, explosive samples or samples capable of forming explosive mixtures.
- ◆ When working with toxic or pathologic substances close hermetically the tubes and strictly observe the safety regulations corresponding to such substances.
- ◆ When you are not using the centrifuge for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the centrifuge by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the centrifuge (see Maintenance).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

6. MAINTENANCE AND CLEANING

To get the best results and a higher duration of this equipment it is essential to follow the processes of use.

Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Balance the centrifuge by using tubes with the same weight, shape and size and put them in geometrically symmetrical positions. If it is necessary, use tubes filled with water. In a centrifuge it is very important the concept of weight so it is advisable to have a balance to compensate the weight of the different tubes.
- ◆ This is absolutely necessary because if the rotor is not well balanced this may produce vibrations that debilitate the centrifuge and adapters increasing the rate of broken tubes and producing a worse sedimentation.
- ◆ Use the appropriate tubes resistant to the R.C.F. used.

Speed adjustment

- ◆ The Nahita Medibas 2740 centrifuge presents an adjustable speed range from 100-4000 r.p.m, in steps of 100 r.p.m. To select the desired speed proceed as the following:
- ◆ Press ▲ ▼ (2.2) buttons, on display the speed value will increase or decrease in steps of 100 r.p.m.

Note: "Speed" display will show at the beginning the speed value selected during last operation.

Time adjustment

- ◆ The Nahita Medibas 2740 centrifuge presents an adjustable time range from 1-99 minutes, in steps of 1 minute. To select the desired time proceed as the following:
- ◆ Press ▲ ▼ (2.4) buttons, on display the time value will increase or decrease in steps of 1 minute.

Note: To work on continuous mode select a 00 time value.

Note: Do not start the centrifuge without closing the lid.

- ◆ Press RUN, the centrifuge starts working.
- ◆ On speed display (2.1), the value of speed will increase until reaching the selected value.
- ◆ On time display (2.3), once the rotor spins at the selected r.p.m., time will decrease (the digit blinks) until reaching 00. If you have selected the continuous mode, time display will show 00 all the time (the digit blinks).

Note: When the centrifuge starts working, the lid is blocked by an electromagnetic lock. Never force its opening with the rotor running.

- ◆ Stop: when the programmed time finishes the rotor will decrease its speed until it stops. When a "click" sounds it means that the lid is unlocked, wait 2-3 seconds before opening making sure the rotor has completely stopped.

Note: When working on continuous mode or if you need to stop the centrifuge before the programmed time, press Stop button.

- ◆ Never try to open the centrifuge before the rotor stops, the lid is blocked.
- ◆ Switch the equipment off by pressing the general/security switch (1.6), to avoid the equipment to start operation accidentally.
- ◆ Now you can take tubes out.

3. TECHNICAL SPECIFICATIONS

Capacity	4 tubes of 100 mL (100x40 mm). 8 tubes Falcon type of 50 mL 16 tubes Falcon type of 15 mL
Speed	1000-4000 r.p.m (adjustable)
F.C.R. max.	3390 g
Timer	1-99 min.
Rotor	Swing-out
Head	Oscillating
Speed regulation	Digital
Time regulation	Digital
Consumption at rest	22 W
Consumption at max r.p.m.	220-400 W
Motor	Elastic suspension with brushes
Environmental conditions of use	5-40° C / RH < 80%
Weigh	28 kg
Dimensions	38x39x37 cm
Fuse	10 A
Supply	220-240 V ± 50-60 HZ

4. THEORETICAL FOUNDATIONS

Centrifugation is a separation technique based on the movement of the particles that are displaced to the distal end of the rotation axis depending on their shape and mass.

The Relative Centrifugal Force (R.C.F.) is the required force to separate particles. It is given as some value times that of gravity (g).


The R.C:F. is calculated with the following formula:

$$F.C.R.= 1.118 \cdot 10^{-5} \cdot r \cdot n^2$$

1,118 * 10⁻⁵ is a constant.

r= radius, the horizontal distance in centimeters from the center of rotation to the bottom of the sample tube.

n= rotation speed expressed in revolutions per minute (r.p.m.).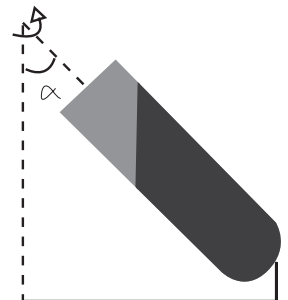

5. INSTALLATION/SETTING UP

Preliminary inspection

Unwrap the centrifuge, take off the involving plastic and take off the polyspan protection in which it comes fitted. Make sure that the equipment does not present any damage because of the shipment. In case it presents any damage, tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.

Note: Since it is very heavy, take the necessary precautions to assure the equipment is properly handled.

Without connecting the centrifuge to the power supply, take out the accessories that are provided inside:

- ◆ Open the lid.
- ◆ With the help of the spanner supplied, loosen the central nut and take out the rotor.
- ◆ Take out the polyspan base with all the accessories.
- ◆ Put again the rotor, by fitting the grooves of the rotor into the bolt of the rotor axe.
- ◆ Put again the washer and the nut and tighten with the spanner.
- ◆ Make sure the rotor is well held and check manually that it turns freely.

Note: During operation, the rotor turns on the opposite direction as the turning direction of the nut; this is to assure that the nut is perfectly tightened and there will be no risk of loosening due to operation.

Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.

Please check that all the accessories are enclosed with the equipment:

- 4 adapters for tubes of 100 ml.
- 4 adapters for tubes of 50 ml.
- 4 adapters for tubes of 15 ml.
- 4 plastic tubes of 100 ml (aprox.)
- 8 falcon standard-like tubes of 50 ml.
- 16 falcon standard-like tubes of 15 ml.
- Spanner.
- Schuko wire.
- Instructions manual.

We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping.

Installation

Before using this instrument, it is convenient for you to familiarize with its components and basic essentials, as well as with its control functions.

PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT WITH THE AIM OF ACHIEVING THE HIGHEST FEATURES AND THE MAXIMUM DURABILITY OF THE INSTRUMENT.


- ◆ Please put the centrifuge on top of a horizontal, plane and stable table making a free space of least at 30 cm per side.
- ◆ Do not put the centrifuge near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc.
- ◆ During operation, people and dangerous material such as flammable or infectious liquids must be out of the working area.
- ◆ The centrifuge is supplied with a standard Shuko wire. Please insert the wire that feeds the AC electric current in the base of current 220V 50Hz $\pm 10\%$ provided with earth wire and to the other end to the centrifuge connector.

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220V 50Hz $\pm 10\%$.

If you are not using the centrifuge for a long period of time please make sure it is disconnected from the net and protected from dust (thus you will avoid accidents and will extend its working-life).

Setting up

- ◆ Open the lid.
- ◆ Place the desire adapters and introduce the tubes inside the buckets. Make sure they are well placed and check that the adapters swing freely. Placing the adapters not properly could produce accidents and damages to the equipment.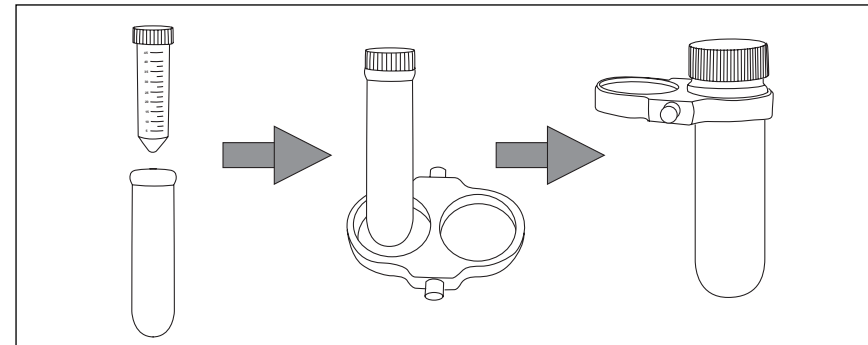

Note: Balance the centrifuge correctly by putting tubes of the same weight, shape and size in position geometrically symmetric; if necessary use tubes filled with water.

- ◆ Close the lid.
- ◆ Switch the centrifuge on by pressing the general/security switch (1.6).