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Discharging liquid

- ◆ Press discharging controller with middle finger, the discharging speed of liquid will be adjusted as per different entering position of filling controller accordingly. When empty the pipette or discharge the liquid with motor is required, please press discharging controller to the end.

**7. FILTER**

The film diameter of this filter is 0.45 µm.

**8. RECHARGEABLE BATTERY**

- ◆ Open the battery cover, push back the knob on the cover with finger, then pull it up.
- ◆ Take out the rechargeable battery.
- ◆ Replacing the new rechargeable battery.
- ◆ Put on the cover, press down the knob and recover.

**9. STORAGE**

- ◆ The stand enclosed can be stuck on the wall and the instrument can be put on it.
- ◆ It can also be put on the table directly

**10. LCD DISPLAY**

- ◆ When the Electromotion Pipette controller stops working for 1 minute, the LCD display will be closed automatically.
- ◆ When pressing the filling controller or “+/-” button, the LCD display will turn on automatically.

**11. CHARGE**

- ◆ When charge up, the charger's indicator is in red, and when it turns to green, the charge is finished.

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At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

Thank you for choosing this equipment. We sincerely wish that you enjoy your Digipette electronic pipette. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

**TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.**

Please bear in mind the following:

- ◆ This manual is inseparable from the Digipette electronic pipette, so it should be available for all the users of this equipment.
- ◆ Follow all required applicable laws about safety and health (electrical security devices, inflammable, toxic and/or pathological products, etc.).
- ◆ You should carefully handle the equipment avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the equipment to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the electronic pipette.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Digipette Technical Assistance Department by email to asistencia@auxilab.es.
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the warranty.
- ◆ Neither fuses nor accessories, including their loss, are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Digipette Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

 **ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

**INDEX OF LANGUAGES**

Spanish ..... 2-5  
 English ..... 6-9  
 French ..... 10-13

**INDEX OF CONTENTS**

1. INTRODUCTION ..... 7  
 2. CONTENTS ..... 7  
 3. PREPARATION PRIOR TO USE ..... 7  
 4. ELECTRICAL POWER SOURCE ..... 8  
 5. DESCRIPTION ..... 8  
 6. OPERATING INSTRUCTION ..... 8  
 7. FILTER ..... 9  
 8. BATTERY ..... 9  
 9. STORAGE ..... 9  
 10. LCD DISPLAY ..... 9  
 11. CHARGE ..... 9

**1. INTRODUCTION**

The electro-motion Pipette Controller is the liquid operating tool which is widely used in Medical and Health, Universities and Colleges and Scientific Research Department. It is simple to use reducing the working intensity greatly, and easily obtaining a satisfactory effect of liquid transfer.

**2. CONTENTS**

Please charge the battery fully before using, it will take about 2-3 hours. Once charge up, it can continuously work for 8 hours.

- Please check the following list when open the box:
- Electomotion Pipette Controller
  - Stand
  - Filter 0.45 μm (2 u.)
  - Charger
  - User's manual
  - Warranty


**3. PREPARATION PRIOR TO USE**

Please charge the battery fully before using, it will take about 2-3 hours. Once charge up, it can continuously work for 8 hours.

**4. ELECTRICAL POWER SOURCE**

- Charger: Input 100-240V-50/60 Hz, Output 5.0V/500mA.
- Battery: Rechargeable Battery 2.4V / 800 mA.

**5. DESCRIPTION**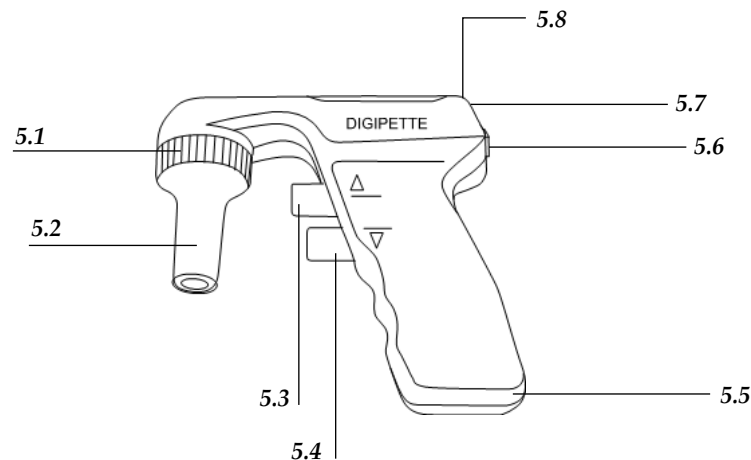

- 5.1 Filter
- 5.2 Silicone cone adapter
- 5.3 Filling controller
- 5.4 Discharging controller
- 5.5 Charge jack
- 5.6 Speed adjusting button
- 5.7 LCD display
- 5.8 Battery

**6. OPERATING INSTRUCTION**

Assembly

- ◆ Insert the pipette
- ◆ Adapt to any size pipette selling in marketing, capacity 1-100 mL

Setting the operating speed

- ◆ Press “+/-” button to adjust the operating speed of the motor as per your demands.
- ◆ When press the filling controller, but the motor no working, press “+” button to increase the motor speed and the working condition will be showed on LCD display.

Filling liquid

- ◆ Press the Filling controller with forefinger, the filling speed of liquid can also be adjusted as per different entering position of filling controller accordingly.