DECLARACIÓN DE CONFORMIDAD CE
CE DECLARATION OF CONFORMITY
DÉCLARATION DE CONFORMITÉ CE

El fabricante | The manufacturer | Le fournisseur:

AUXILAB, S.L.

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◆ Never try to repair the equipment by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...) or damages in nearby equipments.

◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

Cleaning

◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the incubator and produce an early ageing of the equipment.

◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

INSTRUCTIONS ON ENVIRONMENT PROTECTION

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita oven or incubator series 631, 632 or 636. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Please bear in mind the following:

- ◆ This manual is inseparable from the Nahita ovens and incubators, so it should be available for all the users of these equipments.
- ◆ You should carefully handle the oven or incubator avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the oven or incubator to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the oven or incubator.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to asistencia@auxilab.es or by telephone: +34 807 117 040 (0.35 Euros/min).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment's warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

INDEX OF LANGUAGES

Spanish	3-10
English	11-18
French	19-27

INDEX OF CONTENTS

1. USES OF THE INSTRUMENT	12
2. DESCRIPTION	12
3. TECHNICAL SPECIFICATIONS	14
4. INSTALLATION / SETTING UP	15
5. MAINTENANCE AND CLEANING	17
ANNEX I: CE CERTIFICATE	27

1. USES OF THE INSTRUMENT

Nahita ovens and incubators are very useful in research, hospital or industrial laboratories for drying samples, disinfection of material and maintenance of cell and microorganisms cultures.

Made of furnace-painted steel, they present a easy-to-clean stainless steel inner chamber and digital temperature adjustment. To control sample evolution, the series 631 and 632 are provided with a double glass window, while series 636 is equipped with an inner glass door.

2. DESCRIPTION

- 1.1 Control panel
- 1.2 On/off switch
- 1.3 Ventilation knob (only series 631)
- 1.4 Ventilation grilles
- 1.5 Socket

After operation do not forget to turn the equipment off by pressing the on/off switch (1.2).

Security

- ◆ The equipment must be used by previously qualified staff that knows how to operate it thanks to the user manual.
- ◆ Put the equipment on top of a horizontal, plane and stable table making a free space around so as the ventilation grilles are completely free of obstructions.
- ◆ Do not put the equipment near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. Avoid the presence of flammable or toxic products in the working area.
- ◆ During operation dangerous materials such as flammable or pathological substances must be out of the working area.
- ◆ When you are not using the equipment for a long period of time please make sure it is unplugged from the net in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the equipment by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the equipment (see Maintenance).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

5. MAINTENANCE AND CLEANING

To get the best results and a higher duration of the equipment it is essential to follow the processes of use.

Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ◆ The equipment is supplied with a standard Schuko wire. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.

Note: Before the first use or if it has been moved to other location, do not turn the incubator on until 2 hours have passed.

- ◆ Press the on/off switch (1.2).
- ◆ The PV display will show the real temperature inside the equipment, while the SP display will show the previously selected temperature.

Working temperature adjustment

The maximum working temperature is 250 °C for ovens (series 631 and 632) and 80 °C for incubators (series 636). Proceed as follows to select the working temperature:

- ◆ Press Func (2.1). "Sp" will be shown on PV display, indicating that the temperature value can be adjusted.
- ◆ Use the buttons ▲ (2.2) and ▼ (2.3) to select the desired temperature value (shown on SP display). Use the button ◀ (2.4) to go from one digit to the next.
- ◆ Press Func (2.1) twice. The heating resistance will start working until reaching the selected temperature; the real temperature inside the equipment will be shown on PV display, while the SP display will show the selected temperature.

Note: adjust the opening of the ventilation grilles (1.4), manually in series 632 and 636 or by means of the ventilation knob (1.3) in series 631, to adequate it to your needs.


- ◆ If an alarm sounds continuously during operation, it means that the real temperature exceeds the selected temperature in 10 °C for ovens (series 631 and 632) or in 5 °C for incubators (series 636).
- ◆ Let the equipment cool down and set it again.

Working time adjustment


By default, whenever turning the equipment on the working time is off, so the equipment will operate until turning it off by pressing on/off switch (1.2). To work during concrete periods of time proceed as follows:

- ◆ Wait until the equipment reaches the selected working temperature and press Func (2.1) twice. "Sr" will be shown on PV display indicating that the working time can be adjusted.
- ◆ Use the buttons ▲ (2.2) and ▼ (2.3) to select the working time within a range 1-9999 min (shown on SP display). Use the button ◀ (2.4) to go from one digit to the next.
- ◆ Press Func (2.1); the counting down will begin and the SP display will show the remaining working time.
- ◆ Once the selected time has passed, an alarm will sound and Stop will be displayed on SP display. Press Func (2.1) to silence the alarm.

Note: If a new working time has to be programmed, first set time to 0 and press Func (2.1); then select the new working time.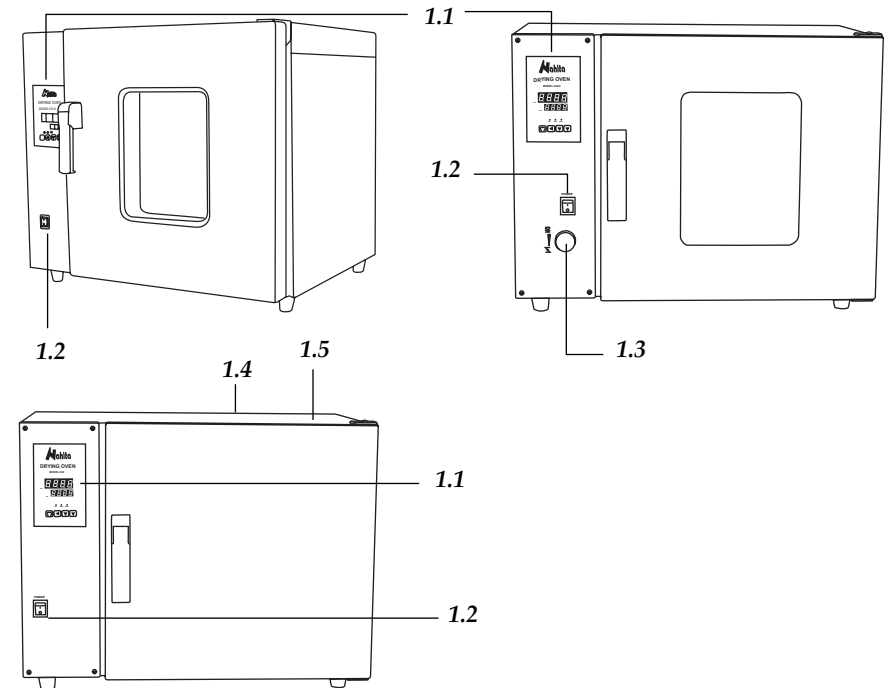

- | | |
|---|---------------------------------------|
| 2.1 Button Func | 2.7 AT indicator (out of use) |
| 2.2 Button ▲ : parameter value increase | 2.8 OUT indicator: resistance working |
| 2.3 Button ▼ : parameter value decrease | 2.9 ALM indicator: alarm |
| 2.4 Button ◀ : go to the next digit | |
| 2.5 PV display | |
| 2.6 SP display | |

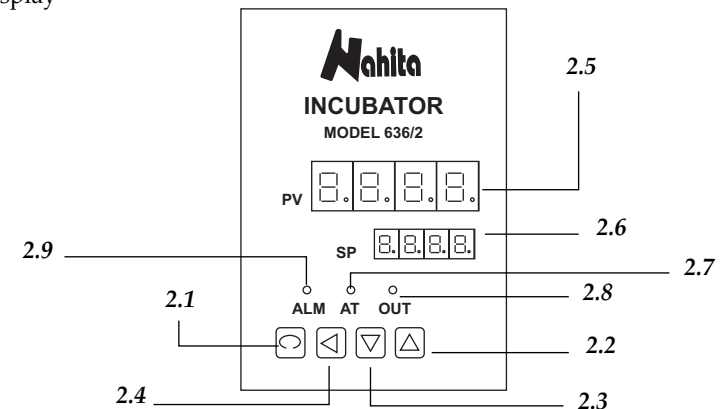3. TECHNICAL SPECIFICATIONS

<i>Series 631 Natural convection</i>	631/2	631/4	631/7	631/13
<i>Code</i>	50631020	50631040	50631060	50631130
<i>Series 632 Forced air</i>	632/2	632/4	632/7	632/13
<i>Code</i>	50632020	50632040	50632060	50632130
<i>Capacity</i>	20 L	40 L	70 L	130 L
<i>Max. temperature</i>	250 °C			
<i>Temperature fluctuation</i>	± 1 °C			
<i>Homogeneity Series 631</i>	3%			
<i>Homogeneity Series 632</i>	2%			
<i>Consumption</i>	1.2 Kw	1.5 Kw	2 Kw	4 Kw
<i>Dimensions (cm)</i>				
<i>External</i>	62x45x52	72x52x59	82x56x69	92x62x79
<i>Internal</i>	28x28x28	35x35x35	45x35x45	55x45x55
<i>Weight</i>	30 Kg	45 Kg	60 Kg	80 Kg
<i>Fuse</i>	15 A			
<i>Power</i>	220 V, 50 Hz ± 10%			

<i>Series</i>	636/2	636/4	636/7	636/13
<i>Code</i>	50636020	50636040	50636060	50636130
<i>Capacity</i>	20 L	40 L	70 L	130 L
<i>Max. temperature</i>	80 °C			
<i>Temperature fluctuation</i>	± 1 °C			
<i>Dimensions (cm)</i>				
<i>External</i>	62x40x52	72x53x57	82x53x68	92x63x78
<i>Internal</i>	28x28x28	35x35x35	45x35x45	55x45x55
<i>Weight</i>	30 Kg	45 Kg	60 Kg	80 Kg
<i>Fuse</i>	15 A			
<i>Power</i>	220 V, 50 Hz ± 10%			

4. INSTALLATION / SETTING UP

Preliminary inspection

◆ Unwrap the equipment, take off the involving plastic and take off the polyspan protection in which it comes fitted. Take off all the protective items and, without connecting the equipment to the net, make sure that it does not present any damage because of the shipment. In case the equipment presents any damage tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.

◆ Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.

◆ Please check that all the accessories are enclosed with the equipment:

- 2 Trays and 4 guides
- Shuko wire
- Warranty
- User's manual

We will only accept the equipment return within 15 days after delivery provided it comes in its original wrapping.

Installation

Before using this instrument, it is convenient for you to familiarize with its components and basic essentials.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

◆ Put the equipment on top of a horizontal, plane and stable table making a free space around so as the ventilation grilles are completely free of obstructions.

◆ Do not put the equipment near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. Avoid the presence of flammable or toxic products in the working area.

◆ Place the trays supplied with the equipment at the desired height.

◆ The equipment is supplied with a Schuko standard wire. Please insert the wire that feeds the AC electric current in the base of current 220 V, 50 Hz ±10% provided with earth wire and the other end to the equipment connector (1.5). The tension should be 220 V, 50 Hz ± 10%.

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220 V, 50 Hz ±10%.