### Agitador basculante | Oscillating stirrer

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Marca|**Trademark:** Nahita

This stirrer is especially suitable for use in laboratories of Haematology thanks to its platform provided with a rocking motion up and down. The stirring speed is soft and constant ensuring an optimal shaking of samples. The platform is coated with a removable rubber tray that avoids tubes to slide.

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*This manual should be available for all users of these equipments. To get the best results and a higher duration of this equipment it is advisable to read carefully this manual and follow the processes of use.*

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Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita oscillating stirrer. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

**TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.**

Please bear in mind the following:

- ◆ This manual is inseparable from the Nahita oscillating stirrer, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the equipment avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the equipment to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the oscillating stirrer.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to [asistencia@auxilab.es](mailto:asistencia@auxilab.es) or by telephone: +34 807 117 040 (0.30 Euros/min).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

ageing of the equipment.

- ◆ The rubber tray can be removed to facilitate cleaning the equipment in case of liquid spillages.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

**ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

#### INSTRUCTIONS ON ENVIRONMENT PROTECTION

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

**Twister  
Medical**

**[www.twistermedical.com](http://www.twistermedical.com)**

**[info@twistermedical.com](mailto:info@twistermedical.com)**

**902 876 983**

- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the stirrer by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the stirrer (see Maintenance).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

## 5. MAINTENANCE AND CLEANING

To get the best results and a higher duration of this equipment it is essential to follow the processes of use.

*Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.*

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ The brushless motor is maintenance-free.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ◆ The stirrer is supplied with a Schuko standard wire. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ◆ Change of fuse: unplugged the stirrer. The fuse compartment is located at the base of the socket. With the help of a sharp object remove the fuse compartment and change the fuse.
- ◆ Never try to repair the stirrer by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the stirrer (burns, hurts...) or damages in nearby equipments.
- ◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

### Cleaning

- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the stirrer and produce an early

**ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

## INDEX OF LANGUAGES

Spanish .....	2-7
English .....	8-13
French.....	14-20

## INDEX OF CONTENTS

1. USES OF THE INSTRUMENT .....	9
2. DESCRIPTION .....	9
3. TECHNICAL SPECIFICATIONS .....	10
4. INSTALLATION / SETTING UP .....	10
5. MAINTENANCE AND CLEANING .....	12
ANNEX I: CE CERTIFICATE .....	20


## 1. USES OF THE INSTRUMENT

This oscillating stirrer with constant speed can be used for several techniques in clinical laboratories of haematology, immunology, etc. and in research laboratories, as for example molecular biology. Thanks to its platform provided with a rocking motion up and down, this instrument assures a soft and uniform homogenization of samples at a constant speed 15 rpm and for unlimited time.

## 2. DESCRIPTION

- 1.1 Switch on/off
- 1.2 Platform with rubber tray
- 1.3 Socket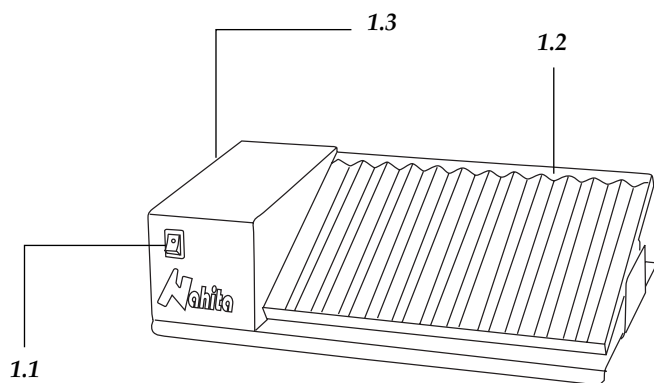

### 3. TECHNICAL SPECIFICATIONS

Code	50683010
Capacity	12 tubes
Max. tube size (ØxL)	20x170 mm
Shaking speed	15 rpm
Platform	255x175 mm
Dimensions (LxWxH)	350x200x95 mm
Weight	2.5 Kg

### 4. INSTALLATION / SETTING UP

#### Preliminary inspection

- ◆ Unwrap the oscillating stirrer, take off the involving plastic and take off the polyspan protection in which it comes fitted. Take off all protections and without connecting the stirrer to the power supply, make sure that it does not present any damage because of the shipment. In case it presents any damage, tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.
- ◆ Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.
- ◆ Please check that all the accessories are enclosed with the equipment:
  - Schuko wire
  - User's manual
  - Warranty certificate

[www.twistermedical.com](http://www.twistermedical.com)

We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping.

#### Installation

Before using the oscillating stirrer, it is convenient for you to familiarize with its components and basic essentials, as well as with its control functions.

**PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT WITH THE AIM OF ACHIEVING THE HIGHEST FEATURES AND THE MAXIMUM DURABILITY OF THE INSTRUMENT.**

- ◆ Put the oscillating stirrer on top of a horizontal, plane and stable table making a free space at least at 30 cm per side. Do not put the equipment near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc.
- ◆ During operation, flammable liquids must be out of the working area.
- ◆ Please insert the Schuko wire that feeds the AC electric current in the base of current 220 V, 50 Hz  $\pm$  10% provided with earth wire and to the other end to the stirrer socket (1.3).

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220 V, 50 Hz  $\pm$  10%.

#### Setting up

- ◆ Put the tubes on the anti-skid grooves perfectly closed to avoid leaks during stirring.
- ◆ Press the on/off switch (1.1) to start with the rocking motion of the stirrer.
- ◆ To stop stirring press again the on/off switch (1.1).
- ◆ If you are not using the stirrer for a long period of time please make sure it is disconnected from the net and protected from dust (this way you will avoid accidents and will extend its working-life).

#### Security

- ◆ The stirrer must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ Put the stirrer on top of a horizontal, plane and stable table making a free space at least at 30 cm per side.
- ◆ Do not put the stirrer near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. Avoid vibrations, dust and dry environments.
- ◆ During operation, flammable liquids must be out of the working area.
- ◆ When you are not using the stirrer for a long period of time please make sure it is unplugged in order to avoid possible accidents.