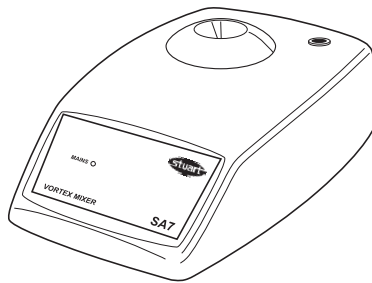
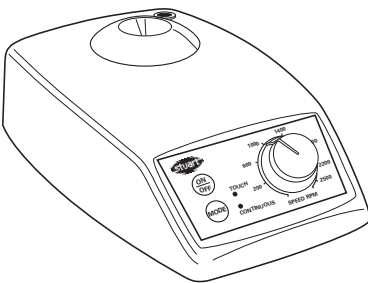
Vortex Mixer SA7 & SA8

Instructions for use
Manuel d'utilisation
Istruzioni per l'uso
Instrucciones de funcionamiento
Bedienungsanweisung

English

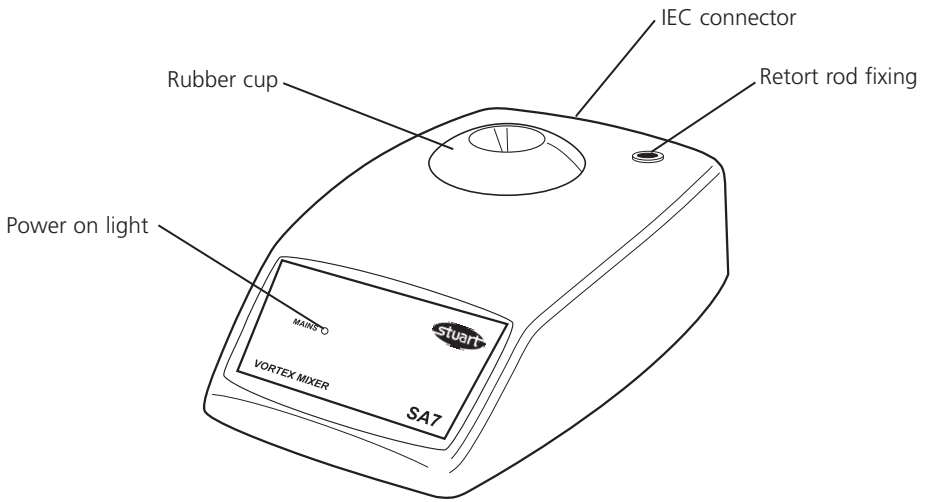
Figure 1: SA7 Fixed Speed model

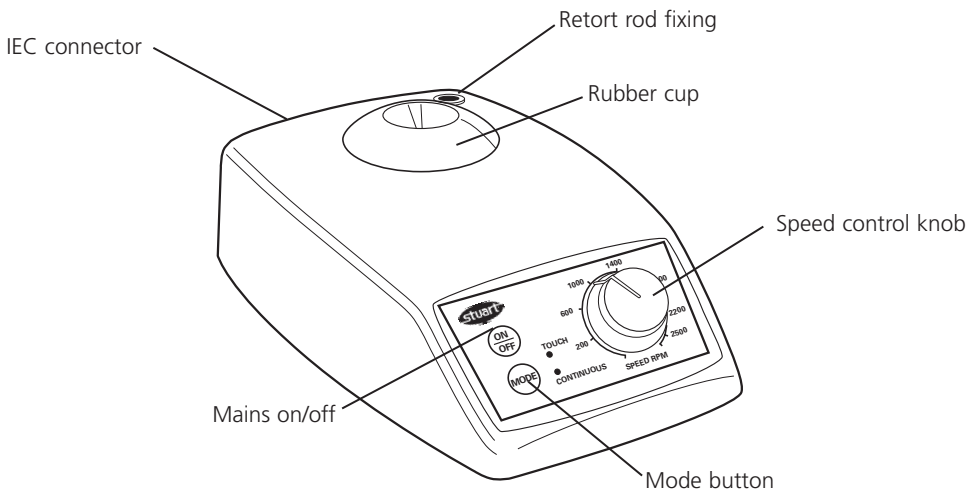
Figure 2: SA8 Variable Speed model

Thank you for purchasing this Bibby Sterilin product. To get the best performance from the equipment, and for your personal safety, please read these instructions carefully before use.

This equipment is designed to operate under the following conditions:-

- ❖ For indoor use only
- ❖ Use in a well ventilated area
- ❖ Ambient temperature range +5°C to +40°C
- ❖ Altitude to 2000m
- ❖ Relative humidity not exceeding 80%
- ❖ Mains supply fluctuations not exceeding 10%
- ❖ Overvoltage category II IEC60364-4-443
- ❖ Pollution degree 2 IEC664
- ❖ Use with a minimum distance all round of 200mm from walls or other items

If the equipment is not used in the manner described in this manual the protection provided by the equipment may be impaired.

Electrical Installation

THIS EQUIPMENT MUST BE EARTHED

Before connection please ensure that the line supply corresponds to that stated on the rating label.

Power requirements:

Model	Wattage
SA7	20W
SA8	20W

There is an IEC socket at the rear of the instrument for connection to the mains supply, (see figure 2). The unit is supplied with two mains leads fitted with IEC plugs for connection to the instrument. One has a U.K. 3 pin plug and the other has a 2 pin "Shuko" plug for connection to the mains. Choose the lead appropriate for your electrical installation and discard the other.

Should neither lead be suitable, take the lead with the U.K. plug and replace the plug with a

suitable alternative. See the enclosed instruction sheet for advice on how to carry out this procedure.

Should the mains lead require replacement a cable of 1mm² of harmonised code H05W-F connected to an IEC 320 plug should be used.

IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN

The mains lead should be connected to the instrument BEFORE connection to the mains supply.

Safety Advice Before Use

- ❖ Never move or carry the unit when in use or connected to the mains electricity supply.
- ❖ Do not mix flammable liquids or use the equipment in hazardous atmospheres.
- ❖ In the case of mains interrupting, a fault or mechanical failure, the unit will NOT continue to operate on removal of fault.
- ❖ Mechanical energy can lead to breakage of glass vessels. Use with care.

Operation

Position the apparatus on a firm level surface, ensuring that the IEC connector is in easy reach.

SA7 Fixed speed version (see Fig 1)

Connect the unit to the mains. Green neon lights. To mix tube, press base firmly down onto the rubber cap, the unit will start. Once the contents of tube are mixed, lift clear, the unit will stop until another tube is pressed onto the cup.

SA8 Variable speed version (see fig 2)

Ensure that the speed control knob is fully to the left. Connect the unit to the mains.

Turn ON by pressing the <On / Off> button once (a green neon lights to show the unit is in the touch mode).

To mix tube, press base firmly down onto the rubber cap. Once the contents of tube are mixed, lift clear. Select the optimum mixing speed by turning the control knob to the right.

To select continuous mode, press the <mode> button (the light will change to show continuous mode). The unit will now remain on until either the <mode> button is pressed to select touch mode or the unit is switched OFF.

Accessory heads

N.B. Accessory heads are for use with model SA8 only. Accessory heads should always be used in the continuous mode.

To change the head, remove the rubber cup by pulling upwards and store in a safe place. Insert the accessory head ensuring that the two slots in the head are in line with the lugs in the vortex mixer top.

The microtitre plate holder (SA8/1) should not be used at speeds above 800rpm.

Maintenance & Servicing

 **WARNING:** Ensure the unit is disconnected from the mains electricity supply before attempting maintenance or servicing.

Periodically clean the instrument using a damp cloth and mild detergent solution. Do not use harsh or abrasive cleaning agents.

Any repairs or replacement of parts MUST be undertaken by suitably qualified personnel.

The following spares and accessories are available from your laboratory dealer.

Description	Catalogue Number
Microtitre plate holder	SA8/1
Spare rubber cup	SA7013
Retort rod, 300mm	SR2

For a comprehensive list of parts required by service engineers conducting internal repairs and a service manual, please contact the Technical Service Department of Bibby Sterilin Ltd. quoting both the model and serial number.

Only spare parts supplied or specified by Bibby Sterilin Ltd. or its agents should be used. Fitting of non-approved parts may affect the performance and safety features designed into the instrument.

If in any doubt, please contact the Technical Service Department of Bibby Sterilin Ltd. or the point of sale.

Bibby Sterilin Ltd
Stone
Staffordshire ST15 0SA
United Kingdom
Tel: +44 (0) 1785 812121
Fax: +44 (0) 1785 813748
E-mail equipment@bibby-sterilin.com
www.bibby-sterilin.com

Warranty

Bibby Sterilin Ltd warrants this instrument to be free from defects in material and workmanship, when used under normal laboratory conditions, for a period of two (2) years. In the event of a justified claim Bibby Sterilin will replace any defective component or replace the unit free of charge.

This warranty does NOT apply if damage is caused by fire, accident, misuse, neglect, incorrect adjustment or repair, damage caused by incorrect installation, adaptation, modification, fitting of non approved parts or repair by unauthorised personnel.

Technical Specification

	SA7	SA8
Speed, rpm	2,500	250-
2,500		
Orbit, mm	4.2	4.2
Continuous mode	No	Yes
Accessory heads	No	Yes
Maximum load		1 Kg
Overall dimension	78 x 135 x 215mm	
Net weight	3.2 kg	
Electrical supply	90 - 240V, 20W	

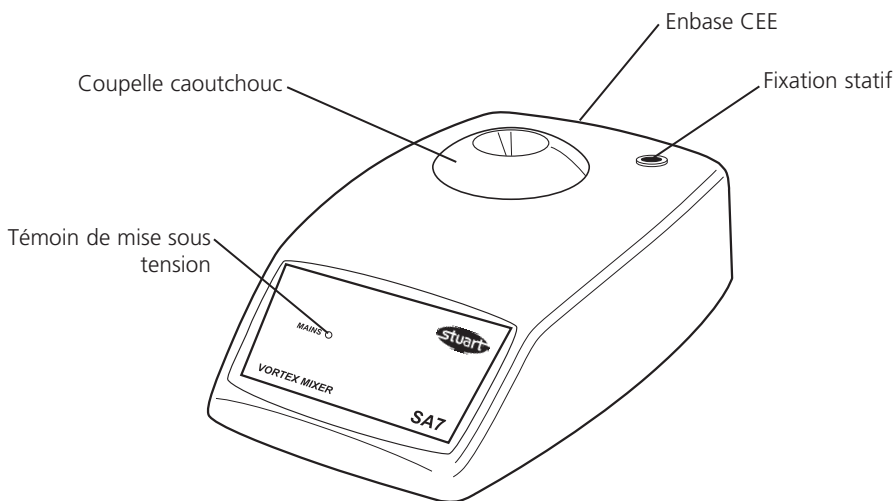
Fig. 1: Version vitesse constante

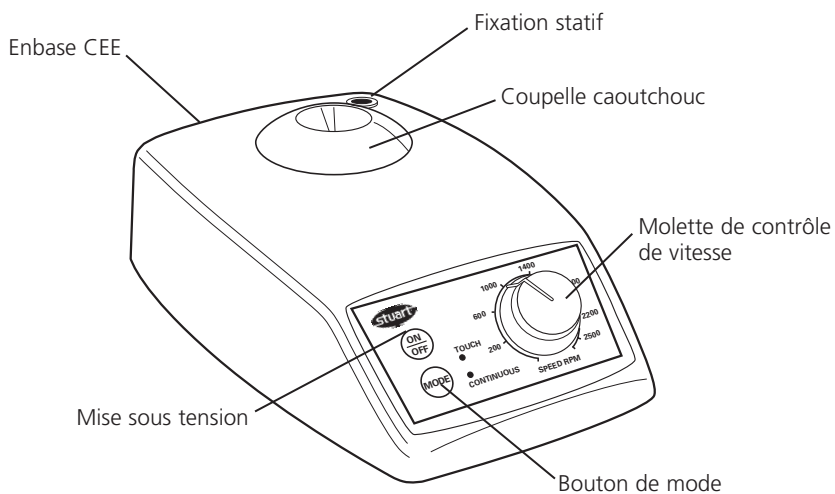
Fig. 2: Version vitesse variable

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Notes

These products meet the relevant EC harmonised standards for radio frequency interference and may be expected not to interfere with, or be affected by, other equipment with similar qualifications. We cannot be sure that other equipment used in their vicinity will meet these standards

and we cannot guarantee that interference will not occur in practice. Where there is a possibility that injury, damage or loss might occur if equipment malfunctions due to radio frequency interference, or for general advice before use, please contact the Technical Service Department of Bibby Sterilin Ltd.

Declaration of Conformity

Model SA7, SA8 VORTEX MIXERS

These products comply with the requirements of the EU Directives listed below:

89/336/EEC Electromagnetic Compatibility Directive, amended by 92/31/EEC and 93/68/EEC.

73/23/EEC Low Voltage Directive amended by 93/68/EEC

Compliance with the requirements of these Directives is claimed by meeting the following standards:

EN61326: 1997, EN61000-4-2: 1995, EN61000-4-4: 1995, EN61000-4-5: 1995
EN61000-4-6: 1996, EN61000-4-11: 1994, EN61000-3-2: 1995, EN61000-3-3: 1995
EN61010-1: 2001, EN61010-2-051: 1996.

The technical documentation required to demonstrate that the products meet the requirements of these Standards is available for inspection on request.

Signed:  (Mr C. R. Baggaley)

Date: 15/8/01

Authority: Technical Manager

for

Bibby Sterilin Ltd - Stone - Staffs - ST15 OSA - UK
Tel: +44 (0) 1785 812121 - Fax +44 (0) 1785 813748

INSPECTION REPORT

MODEL _____

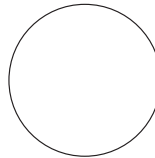
ELECTRICAL SAFETY

- 1. Earth continuity
- 2. Insulation
- 3. Flash test

FUNCTIONAL

- 1. Indicators
- 2. Vortex action
- 3. Visual acceptance

QUALITY CONTROL INSPECTOR

Bibby Sterilin

Bibby Sterilin Ltd

Tilling Drive Stone
Staffordshire ST15 0SA
United Kingdom
Tel: +44 (0)1785 812121
Fax: +44 (0)1785 813748
www.bibby-sterilin.com

Bibby Sterilin France

BP 79-77793 Nemours Cedex
France
Tel: +33 1 64 45 13 13
Fax: +33 1 64 45 13 00
bibby@bibby-sterilin.fr

Bibby Sterilin Italia

Via Alcide de Gasperi 56
20077 Riozzo Di Cerro Al Lambro
Milano
Italy
Tel: +392 983 8969
Fax: +392 982 30211

Afora S.A.

Calle Aribau 240
08006 Barcelona
Spain
Tel: +343 93-306 98 00
Fax: +343 93-306 98 23
e-mail: marketing@afora.com
www.afora.com