**BAÑO ULTRASONIDOS
ULTRASONIC WATER BATH
BAIN ULTRASONS**

Nahita

Modelo / Model / Modèle 620-1

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5. MAINTENANCE ET NETTOYAGE14

Security

- ◆ The ultrasonic bath has to be used by qualified personal so it gets its correct use.
- ◆ You should put the bath in a horizontal plane stable table, having a free space of at least 30 cm per side.
- ◆ Do not put the bath near any warm supply (burners, blowlamp...), do not expose directly to the sun, avoid vibrations and dust. During its functioning, do not use the equipment near fluids that can give off steams or produce explosive or inflammable mixtures.
- ◆ If you are not going to use the equipment during a long period of time please make sure that it is unplugged from the net.
- ◆ Never dismantle the different pieces of the ultrasonic water bath to repair yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty.
- ◆ Fabricated under European regulations.

5. MAINTENANCE & CLEANING

To get the best results and a higher duration of this equipment it is necessary to follow the processes of use.

Note: All the processes of use mentioned before will not have any value unless you keep a continued and careful maintenance.

- ◆ Please follow the processes of use of this manual
- ◆ This manual should be available for all users of these equipments
- ◆ Always use original components and supplies. Other dispositives can be similar but they can damage the equipment.
- ◆ The bath is suminstrated with a Schuko standard wire and it should be connected to a current wire provided with an earth wire. It should be handy to be disconnected in case of emergency.
- ◆ In case of damage or doubt it can be solved getting into touch with the NAHITA Technical Service
- ◆ For any manipulation or cleaning of the equipment please make sure it is disconnected from the net.
- ◆ Never dismantle the different pieces of the ultrasonic water bath to repair yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty.

Cleaning

- ◆ For the cleaning of the inox steam parts use always a special cloth that does not give off fluff, do not use any cloth that can produce damages to the equipment.
- ◆ For the cleaning of the different pieces of the equipment , we highly recommend to use neutral products that do not have a PH superior than 6 (soaped water without abrasives).

ATTENTION!! IF AN EQUIPMENT IS NOT PROPERLY CLEAN AND DESINFECTED IT WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE

Thank you for choosing this equipment. We sincerely wish that you enjoy your Ultrasonic Water Bath NAHITA, we highly recommend to look after this equipment according to what is stated in this manual.

NAHITA develops its products being remarkable in the ergonomoy and security of its user. The incorrect use of the equipment, can cause accidents and electric discharges, circuit breakers, fires, damages, please read the point of maintenance , where we expose the security notes.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

TO GET THE BEST RESULTS AND HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Please bear in mind the following:

- ◆ This manual should be available for all the users of these equipments
- ◆ Please use always the plastic baskets placing the articles you are going to clean, avoiding them to contact with the stainless steel tank, this way we will preserve the life of the equipment.
- ◆ We should manipulate carefully avoiding the sudden movements, knocks, free fall of heavy objects, manipulation of sharp objects.
- ◆ Never dismantle the different pieces of the ultrasonic water bath to repair yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty.
- ◆ Do not use the equipment with fluids that can give off steams or produce explosive or inflammable mixtures.
- ◆ If you have any doubt do not hesitate in contacting your distributor or our NAHITA Technical Assistant Department by email asistencia@auxilab.es, by telephone 34 948310513 or by fax 34 948310500 or 948312071.
- ◆ Please do not forget filling the Warranty Certificate and send it before 15 days after the date of purchase.

ATTENTION!! IF AN EQUIPMENT IS NOT PROPERLY CLEAN AND DESINFECTED IT WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE

INDEX OF CONTENTS

1. USES OF THE INSTRUMENT	8
2. DESCRIPTION	8
3. TECHNICAL SPECIFICATIONS.....	8
4. SETTING UP	9
5. MAINTENANCE & CLEANING.....	10


1. USES OF THE INSTRUMENT

The cleaning by Ultrasonic baths consists in the use of the sound waves of high frequency (40/50Hz) they make in the interior of the tin. A cavitation effect is produced and causes the formation of millions of microscopic bubbles of low pressure that make a molecular cleaning, eliminating the impurities, pollutants and dirt of the pieces or materials that must be cleaned.

The Ultrasonic Nahita cleaning baths allow all kind of instrument cleaning without detaching them. Soffidicated components or those that have a special difficulty to be cleaned, can be cleaned without being separated, thanks to the cavitation going in to every item where the cleaner solution touches the surface.

They are suitable for: odontology and surgical material, as well as in jewelry, clockwork, optics and small tools, all kind of plastic, glass, porcelain, metal materials...

2. DESCRIPTION

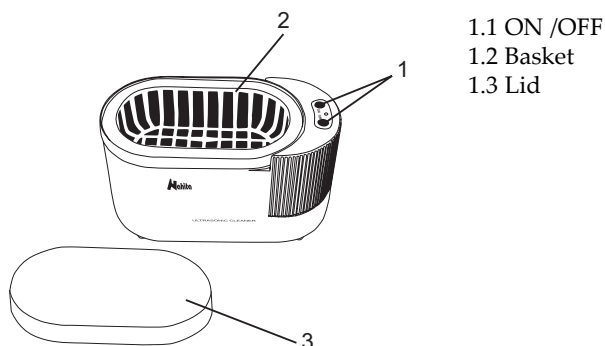3. TECHNICAL SPECIFICATIONS

Feeding	220 V \pm 10%, 50-60Hz
Frequency	40KHz
Dimensions	223 x 114 x 123 mm
Inner size	167 x 90 x 52 mm
Volume	0.6 l
Weight	1Kg

The ultrasonic bath is made of special plastic and a stainless steel tank, it also has a plastic basket where you can deposit the pieces for their cleaning, it also has a plastic lid to avoid the splashing.

4. SETTING UP

Preliminary inspection

- ◆ Unrrape the bath, take off the envolving plastic and take off the poliespan protection in wich it comes fitted. Please make sure that it does not present any damage because of the shipment. If it does please communicate to your transport agent or suministrator immediately, so that they can make the claims in the correct time limit.
- ◆ Please keep the original wrapping, you will always need it for returns enclosed with all the accesorize suministrated.
- ◆ Please check that all the accesorize are enclosed with the equipment:
 - Plastic basket
 - Plastic lid
 - Instruction Manual
 - Warranty Certificate

Installation

- ◆ Before using this instrument, it is convinient for you to familiarize with its components and basic essentials.

TO GET THE BEST RESULTS AND HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

- ◆ Please put the bath on top of a horizontal, plane, stable table making a free space at least at 30 cm per side of the bath. Do not put the bath near any warm supply (burners, blowlamp...), do not expose directly to the sun, etc. During its functioning, do not use the equipment near fluids that can give off steams or produce explosive or inflammable mixtures.
- ◆ The bath is suministrated with a Schuko standard wire.
- ◆ Please insert the wire that feeds the A/C in the base of current 220V \pm 10% 50-60Hz provided with earth wire.

The manufacturer and the distributor will not assume any responsibility for the damages produced to the equipment in its installing or damages to persons suffered by the wrong use of the electric connection. The tension should be 220V \pm 10% 50-60Hz.

- ◆ If you are not using the bath during a long period of time please make sure it is disconnected from the net.

Setting up

- ◆ The bath has a Schuko cable for the net; it should be connected to a current wire provided with an earth wire. It should be handy to be disconnected in case of emergency.
- ◆ For covering the objects you want to clean, you have to introduce a sufficient quantity of water. You also have to add a soap solution for the cleaning with the ultrasonic (free from phosphates), then you proceed to connect to the current for its use.
- ◆ Turn on the equipment pressing the button ON. Once enough time has passed to clean the object turn off the equipment by the button OFF. If the time programed by the clock has passed it will turn off automatically (this time is around 15 minutes). You can repeat the times you need this cycle for the correct cleaning of the objective. You can also interrupt it (OFF) in case you need less time than programmed.