**AGITADOR MAGNÉTICO DISPLAY LED 20 L
DISPLAY LED MAGNETIC STIRRER 20 L
AGITATEUR MAGNETIQUE DISPLAY LED 20 L**

Modelo / Model / Modèle / RSLAB-3C

RS Lab

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**INSTRUCTIONS ON ENVIRONMENT PROTECTION**

At the end of its life cycle, please, does not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non-adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

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INDEX OF CONTENTS

1. Safety instructions	10
2. Proper use	11
3. Inspection	11
4. Trial run	11
5. Control and display	12
6. Faults	13
7. Maintenance and Cleaning	13
8. Associated Standards and regulation	14
9. Technical data	14
10. Accessories	15

Welcome to the magnetic stirrer RSLAB-3C (heating) user manual. Users should read this manual carefully, follow the instructions and procedures, and beware of all the cautions when using this instrument.

When help needed, you can always contact the service department of manufacturer or your supplier for technical support. Please provide the customer care representative with the following information: Serial number, description of problem, methods and procedures adopted to resolve the problems and your contact information.

These RSLAB stirrers are warranted to be free from defects in materials and workmanship under normal use and service for a period of 24 months from the date of invoice. The warranty is extended only to the original purchaser. It shall not apply to any product or parts which have been damaged on account of improper installation, improper connections, misuse, accident or abnormal conditions of operation.

For claims under the warranty please contact your local dealer. You may also send the instrument direct to our works, enclosing the invoice copy and giving reasons for the claim.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

1- SAFETY INSTRUCTIONS

Risk of burn!

♦ Caution when touching the housing parts and the heating plate. The heating plate can reach temperatures of 550 °C.

♦ Pay attention to the residual heat after switching off.

Protective ground contact!

♦ Make sure that socket is earthed (protective ground contact) before use.

♦ When work, wear the personal guard to avoid the risk from splashing and evaporation of liquids and release of toxic or combustible gases.

♦ Set up the instrument in a spacious area on a stable, clean, non-slip, dry and fireproof surface, do not operate the instrument in explosive atmospheres, with hazardous substances or under water.

♦ Gradually increase the speed, reduce the speed if the stirring bar breakaway because of too high speed and if the instrument is not running smoothly, or if the container moves on the base plate.

♦ Temperature must always be set to at least 50 °C lower than the fire point of the media used.

♦ Beware of hazards due to flammable materials or media with a low boiling temperature, overfilling of media and unsafe container.

♦ Process pathogenic materials only in closed vessels.

♦ Check the instrument and accessories before handle it for damaged component. Safe operation is only guaranteed with the accessories described in the "Accessories" chapter. Accessories must be securely attached to the device and cannot come off by themselves. Always disconnect the plug before fitting accessories.

♦ Ensure that the external temperature sensor is inserted in the media to a depth of at least 20 mm.

♦ When using metal vessels, do not place the temperature sensors on the bottom of the vessel. Placing sensors on the vessel bottom can cause excessively high temperatures to be measured especially in media which have poor conductivity. The tip of the measuring sensor must be at least 5 mm from the vessel bottom; a distance of 10 mm is ideal.

♦ The instrument can only be disconnected from the main power supply by pulling out the main plug or the connector plug.

♦ The voltage started on the label must correspond to the main power supply.

♦ Ensure that the main power supply cable does not touch the heating base plate. Do not cover the device.

10- ACCESSORIES (NOT INCLUDED)

[1] 90694001

[2] 90694002

[3] 90694011

[4] 90694305

[5] 90694306

[6] 90694307

[7] 90694308

[8] 90694309

[9] 90694310

Reference

- [1] 90694001
- [2] 90694002
- [3] 90694011
- [4] 90694305
- [5] 90694306
- [6] 90694307
- [7] 90694308
- [8] 90694309
- [9] 90694310

Description

- Temperature sensor PT1000, l. 230 mm
- Stand kit for sensor PT1000
- Temperature sensor PT1000, with glass coated, l. 230 mm
- Poly-bloc stand red, 11 holes ø 12,5 mm / width 20 mm / tubes 4 ml
- Poly-bloc stand purple, 4 holes ø 28 mm / width 24 mm / tubos de 20 ml
- Poly-bloc stand blue, 4 holes ø 28 mm / width 30 mm / tubos de 30 ml
- Poly-bloc stand black, 4 holes ø 28 mm / width 43 mm / tubos de 40 ml
- Poly-bloc stand green, 6 holes ø 17,8 mm / width 26 mm / tubos de 8 ml
- Poly-bloc stand gold, 4 holes ø 21,6 mm / prof. 31,7 mm / tubos de 16 ml

8- ASSOCIATED STANDARDS AND REGULATIONS

Construction in accordance with the following safety standards:

- EN 61010-1
- UL 3101-1
- CAN/CSA C22.2 (1010-1)
- EN 61010-2-10

Construction in accordance with the following EMC standards:

- EN 61326-1


Associated EU guidelines:

- EMC guidelines: 89/336/EEG
- Instrument guidelines: 73/023/EEG

9- TECHNICAL DATA

<i>Voltage</i>	100-120/200-240 VAC
<i>Frequency</i>	50/60 Hz
<i>Power</i>	50-1050W (heating model)
<i>Stirring point position quantity</i>	1
<i>Max. stirring quantity (H2O)</i>	20 Liters
<i>Max. magnetic bar</i>	80x10 mm
<i>Max. power input of motor</i>	18 W
<i>Max. power output of motor</i>	10 W
<i>Speed range</i>	0-1500 rpm
<i>Rotary speed display</i>	Scale
<i>Speed display accuracy</i>	1 rpm
<i>Hotplate material</i>	Glass ceramic
<i>Dimensions of workplate</i>	184x184 mm
<i>Dimensions of heating</i>	176x176 mm
<i>Heating power</i>	1000 W
<i>Temperature range</i>	RT-550 °C
<i>Temperature display</i>	LED
<i>The safety temperature range of hot plate</i>	100 a 580°C
<i>Dimensions</i>	215x360x112 mm
<i>Weight</i>	5,3 Kg
<i>Permitted ambient temperature</i>	5-40°C
<i>Permitted relative humidity</i>	80%
<i>Protection class ac. to DIN 60529</i>	IP21
<i>RS232 Interface</i>	No

- ◆ The instrument may only be opened by experts.
- ◆ Keep away from high magnetic field.
- ◆ Observe the minimum distances between the devices, between the instruments and the wall and above the assembly (min. 100 mm).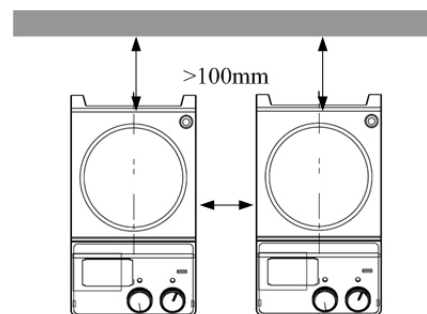

2- PROPER USE

The instrument is designed for mixing and/or heating liquids in schools, laboratories or factories. This device is not suitable for using residential areas or other constraints mentioned in chapter 1.

3- INSPECTION

- ◆ Unpack the equipment carefully and check for any damages which may have arisen during transport. If it happens, please contact manufacturer for technical support.
- ◆ If there is any apparent damage to the system, please do not plug it into the power line.
- ◆ The RSLAB-3C includes the following items:
 - Main unit: 1 unit
 - Power cable: 1 unit
 - Stirrer bar: 1 unit
 - User manual: 1 unit


4- TRIAL RUN

- ◆ Make sure the required operating voltage and power supply voltage match.
- ◆ Ensure the socket must be earthed reliably.
- ◆ Plug in the power cable, ensure the power be on and begin initializing.
Add the medium into the vessel with a stirring bar.
- ◆ Put the vessel on the workplate.

- ◆ Set the rated stirring speed and start stirring.
- ◆ Set the rated temperature and start heating.
- ◆ If these operations above are normal, the device is ready to operate. If these operations are not normal, the device may be damaged during transportation, please contact manufacturer for technical support.

5- CONTROL AND DISPLAY

1. Display
2. LED heat display
3. LED stir display
4. Heating knob
5. Stirring knob
6. Switch ON/OFF
7. Safe Temperature Button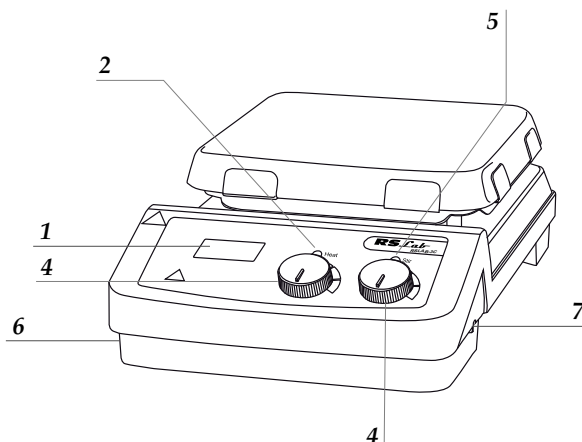

Items	Description
Stirring Knob	Set the rated rotary speed from 0 to 1500 rpm. The stirring function is switched on or off by pushing the knob.
Stirring Knob	Set the rated temperature. The heating function is switched on or off by pushing the knob
LED Heat	When switched on, the LED heat is lit
LED Stir	When switched on, the LED stir is lit
Safe Temperature Button	Set safety temperature, stop heating when exceed secure temperature
Power Switch	Switch on or off the instrument
Display LED	Display of the selected and actual heating T

6- FAULTS

- ◆ Device can't be power ON:
 - Check whether the power cable is plugged
 - Check whether the fuse is broken or loose
- ◆ Fault in power on self-test:
 - Switch OFF the unit, then switch on and reset the instruments to factory default setting
- ◆ Temperature cannot reach set point:
 - Check whether the safety temperature value is set too low
- ◆ Stir speed cannot reach the set point:
 - Check whether the stirring bar breakaway monitoring is on
 - Excessive medium viscosity may cause abnormal speed reduction of the motor
- ◆ Heating or stirring cannot be started when push the temperature control knob:
 - Check whether the unit is in the mode setting program
 - Exit from the mode setting and restart the heating/stirring functions

If these faults are not resolved, please contact manufacturer for technical support.

7- MAINTENANCE AND CLEANING

Proper maintenance can keep instruments working in a good state and lengthen its lifetime.

- ◆ Be careful not spray the cleaners into the instrument when cleaning.
- ◆ Unplug the power line when cleaning.
- ◆ Only use cleaners that we advise as below:
 - Dyes: isopropyl alcohol
 - Construction materials: water containing tenside / Isopropyl alcohol
 - Cosmetics: water containing tenside / Isopropyl alcohol
 - Foodstuffs: water containing tenside
 - Fuels: water containing tenside
- ◆ Wear proper protective gloves during cleaning of the instrument.
- ◆ Before using other method for cleaning or decontamination, the user must ascertain with the manufacturer that this method does no destroy the instrument.
- ◆ Unit should be clean and fitted in the original packing or a suitable packing for transport.
- ◆ The enamel makes the heating plate easier to care for and more resistant to acids and bases. Because of it, however, the heating plate is also more susceptible to extreme fluctuations in temperature and the force of impact. This can result in cracks forming or the coating flaking off.