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Thanks for choosing a NAHITA Electronic Balance, its high mechanical quality will allow you to make use of this equipment for a long years. Nevertheless to get the best results and a higher duration out of this instrument it is highly advisable to read this manual and follow the instructions of use.

BEFORE CONNECTING, OPERATING THIS INSTRUMENT, READ THESE INSTRUCTIONS THOROUGHLY.

In order to use this equipment, bear in mind the following points:

- ◆ This manual is inseparable from the balance and must be available for all users of this equipment.
- ◆ Any doubts can be clarified by getting in touch with the NAHITA service.
- ◆ As an accuracy instrument, it should be handled carefully avoiding the sudden movements and knocks.
- ◆ Never dismantle the different pieces of the balance to repair them yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty.

ATTENTION !! NO EQUIPMENT WILL BE REPAIRED BY OUR TECHNICAL SERVICE UNLESS IT IS PROPERLY CLEAN AND DISINFECTED

INDEX OF CONTENTS

1. APLICACIONES.....	9
2. DESCRIPTION.....	9
3. TECHNICAL SPECIFICATIONS.....	10
4. INSTALATION.....	10
5. FUNCTIONS.....	12
6. MAINTENANCE AND CLEANING.....	12

- ◆ Save all packing material until you are sure that the instrument works correctly. All defective items must be returned in the original packing together with the supplied accessories.
- ◆ It is prohibited to use strong chemical solvent to clean the surface.
- ◆ Do not leave the balance exposed to direct sunlight for long periods of time. This may deform or discolor the cabinet and may also cause malfunctioning.
- ◆ Please take out the dry batteries if it's to be left unused for a long time, wipe the balance, enclosed it with dryer in a plastic pocket.
- ◆ Do not press the display, Do not press the front panel with sharp objects.
- ◆ Don't dip the balance into water, don't leave it out in the rain or even don't wash it with water. In case some water drops, please dry it immediately with a cloth.

INSTRUCTIONS ON ENVIRONMENT PROTECTION

Do not throw out this balance in the ordinary garbage; take it to a collection point for recycling electrical and electronic equipments. Does not contain dangerous or toxic elements to humans.

Materials are recyclable. Recycling is an important contribution to environmental protection. Please contact the administration of your community for be advised on the collection points.

5. DIFFERENT FUNCTIONS OF THE BALANCE

Tare function

- ◆ Tare function is used when you want to weigh any substance without but not the recipient containing it. For instance, when you want to weigh liquids or solid reagents that you are going to dissolve afterwards. You should do the following:
 - ◆ Make sure there is nothing on the pan (1.15) and check the display shows value 0 or 0.0 (depending on gradation you choose).
 - ◆ Put the recipient on the balance's pan (1.15) and then, press TARE key (1.5). After this, you will read 0 or 0.0 on the weight display (depend gradation you choose). The balance will be now ready to start weighing.
 - ◆ Once the weighing process is finished, remove the recipient from the balance. You will be able to read the recipient's weight with a negative value on the display.
 - ◆ To recover its usual mode press TARE key (1.5) again. The display should show 0 or 0.0 again.

Piece counting function

- ◆ Make sure there is nothing on the pan (1.15), check weight display (1.3) reads 0 or 0.0 (depend gradation you choose). If you need to put some recipient you will use tare function.
- ◆ Place a unit of product to be counted. Bear in mind sensitivity that will depend on the balance model. If the weight of the piece is smaller than indicated, then, place 10 units. When the balance is stable, introduce the unit weight by pressing the numbers keys (1.6). On the unit weight display, you can read the weight as you press the key. Press ACCU(1.11), Remove the piece.
 - ◆ The balance is now ready to start counting. Place the product you want to count. On displays you will read:
 - Number display (1.1) shows the number piece.
 - Unit Weight display (1.2) shows the weight piece.
 - Weight display(1.3) shows the total weight piece.
 - ◆ Once finished the work, for cleaning data, press CLEAR key (1.6), press ACCU key(1.8) and them wait for some seconds. Press again CLEAR key(1.6). Now, previous data is deleted and you can start again.

6. MAINTENANCE AND CLEANING

- ◆ Preserve the balance from heavy shocks and look for a dry place, away from dust, direct sun beams or air flows. Handle the balance gently, since it is an accurate instrument. Use smooth gauze with a soapy solution to clean it.
- ◆ Never remove its mechanical components. You will not only lose your warranty, but also you can damage your balance seriously.
- ◆ Don't throw away the package. It might be useful in case you need to move it, to send it if it needs repairing or you stop using it for a long time in the future.


1. INTRODUCTION

NAHITA electronic balances are basic instruments in weighing tasks in every laboratory. They have been developed with the latest technology. They are of a great reliability and durability. Their solid, clean and ergonomic design is very suitable for any industrial usage. It's smooth surface makes the cleaning easier.

They are designed for every kind of uses in lots of fields thanks to the function to count pieces. They are used in teaching, laboratories, pharmacy and industry, chemistry, etc...

2. DESCRIPTION

- | | |
|------------------------|--------------------------------|
| 1. Number Display | 9. SAMP key |
| 2. Unit weight Display | 10. Power light display. |
| 3. Weight Display | 11. Power ON/OFF |
| 4. ZERO key | 12. Numer Key |
| 5. TARA key | 13. Indicator of AC connection |
| 6. CLEAR key | 14. Level |
| 7. SET key | 15. Platform. |
| 8. ACCU key | |

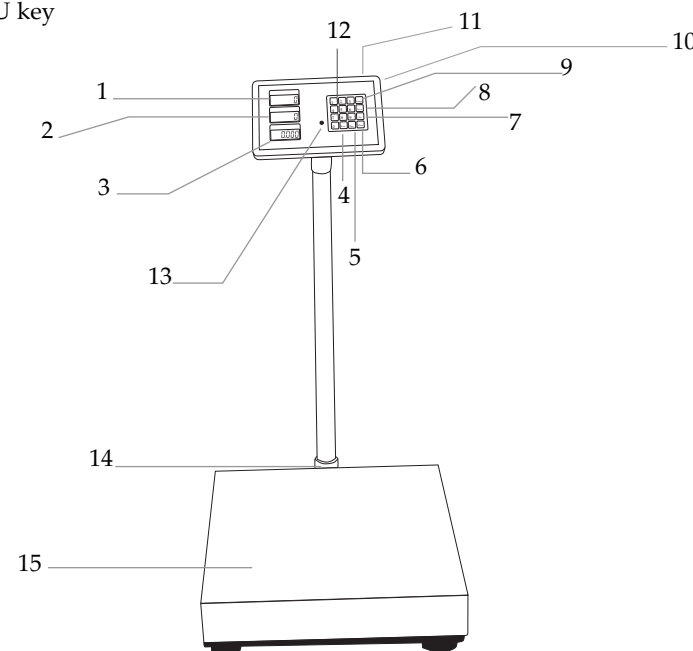3. TECHNICAL SPECIFICATIONS

<i>Code</i>	55059460	55059510	55059530
<i>Capacity</i>	60 Kg	150 Kg	300 Kg
<i>Sensibility</i>	2 g	5 g	10 g
<i>Platform dimensions</i>	450x600 mm.		
<i>Time of stabilization</i>	10 segundos		
<i>Working temperature</i>	5 a 40° C		
<i>Power supply</i>	220V, 50 HZ		

4. OPERATING INSTRUCTIONS

Preliminary examination

- ◆ Remove the instrument from the packing material and examine it carefully to make sure no damage has occurred during shipping. If there is any damage, notify to your dealer.
- ◆ Keep all packing material until you are sure that the instrument works correctly. All defective items must be returned in the original packing together with any other supplied accessories.

Installation

Before using this instrument, please carefully read this manual for its correct use.

BEFORE CONNECTING, OPERATING THIS INSTRUMENT, READ THESE INSTRUCTIONS THOROUGHLY.

- ◆ The instrument must be placed on a flat, horizontal desk. Avoid wetness, heat, vibrations and airflows. The environmental lighting must not be too bright. Keep away from any magnetic force, electric field or gadget with can produce a high frequency field.
- ◆ Place the pan on its place, it must be inserted in the four holes that the balance has for this purpose.
- ◆ Check the level the balance incorporates in its back side is in the appropriate position (in the centre). If it is in a wrong position, try centering the bubble with the knobs close to the balance legs in the front and rear parts.

Insert the tube in the orifice that is in the part down of platform, Subject with the screw. Place the control panel in the last of the tube and subject with the screw. Connect the plug of 9 pin that has the tube, in the control panel.

Charging the batteries

- ◆ Before using the balance, its battery must be completely charged.
- ◆ In order to do so, insert the supplied cable in the schucko pin, the plug it in to an inta-

ke earth connection

- ◆ The battery will be charged after approximately 20 h of connection. Note: the indicator of AC connection (13) will be continuously on while the equipment is connected to the power supply.

Neither the manufacturer nor the distributor will take any responsibility on damaged equipments or personal injuries due to an inappropriate connection to electricity.

Environmental work conditions

<i>Working temperature range</i>	5°-40° C
<i>Relative humidity</i>	50-85%

Starting

- ◆ Switch on the power ON/OFF(1.11), on the right-hand side
- ◆ Check the level the balance incorporates in its back side (1.14) is in the appropriate position (in the centre). If it is in a wrong position, try centering the bubble with the knobs close to the balance legs in the front and rear parts.
- ◆ Balance begin self-testing until getting normal weighting condition. The displays will turn on and the balance will carry out a selftesting.
- ◆ To choose the graduation/sensibility you want to work with use the SET key(1.7). Model capacity divided into graduation gives the weighing sensibility. The higher graduation, the better sensitivity the balance will work with. Look at table.

<i>Code</i>	<i>Capacity</i>	<i>Graduation</i>	<i>Sensibility</i>
55050360	60 g	3000	20 g
		6000	10 g
		12000	5 g
		30000	2 g
55050410	150 g	3000	50 g
		6000	20 g
		12000	10 g
		30000	5 g
55050430	300 g	3000	100 g
		6000	50 g
		12000	20 g
		30000	10 g

- ◆ The balance will be ready to start weighing.
- ◆ Turn off (Press ON/OFF) (1.11)

Try recharging battery once it is completely discharged, this way duration of battery will increase; LB (Low Battery) indicator will be displayed only when the equipment is disconnected from the power supply.