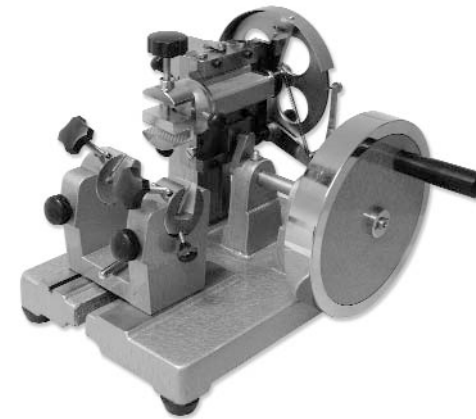
Modelo / Model / Modèle 511

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Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita microtome model 511. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user. The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of *Maintenance*, where we expose the security notes.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Please bear in mind the following:

- ◆ This manual is inseparable from the microtome nahita model 511, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the microtome avoiding sudden movements, knocks, free fall of heavy / sharp objects on it.
- ◆ Never dismantle the different pieces of the microtome to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the microtome.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. you can also tell us any doubts or suggestions you have by contacting nahita technical assistance department by email to asistencia@auxilab.es or by telephone: +34 807 117 040 (0,30 euros/min).
- ◆ This equipment is protected under the warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the microtome warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the microtome warranty.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. in case you have to send the equipment to nahita technical assistance department you should enclose the original invoice or a copy as guarantee.
- ◆ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

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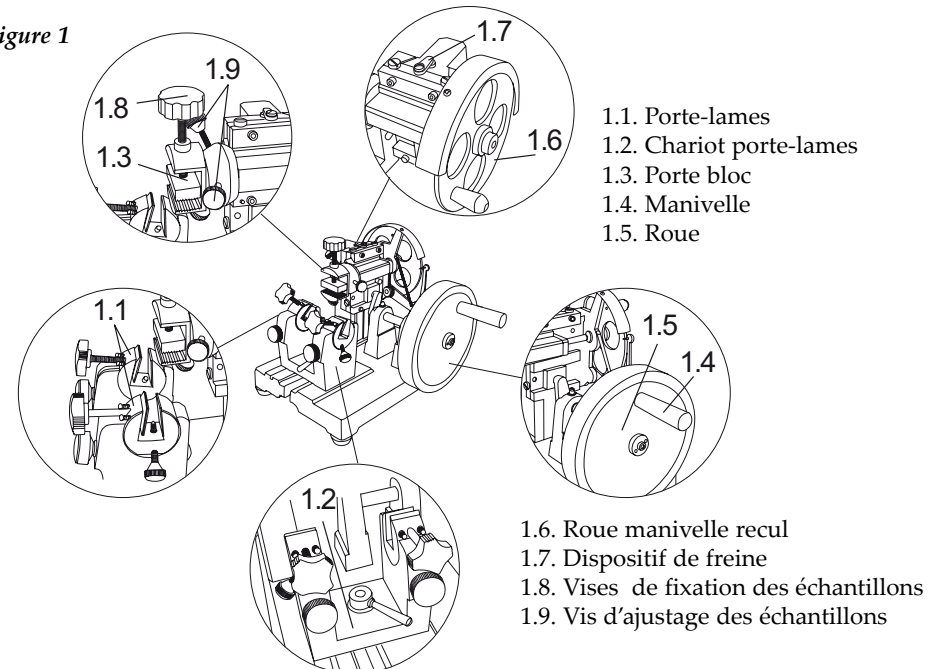
INDEX OF CONTENTS

1. USES OF THE INSTRUMENT	11
2. DESCRIPTION	11
3. TECHNICAL SPECIFICATIONS	12
4. INSTALLATION / SETTING UP	12
5. MAINTENANCE AND CLEANING	16
ANNEX I: CE CERTIFICATE	17

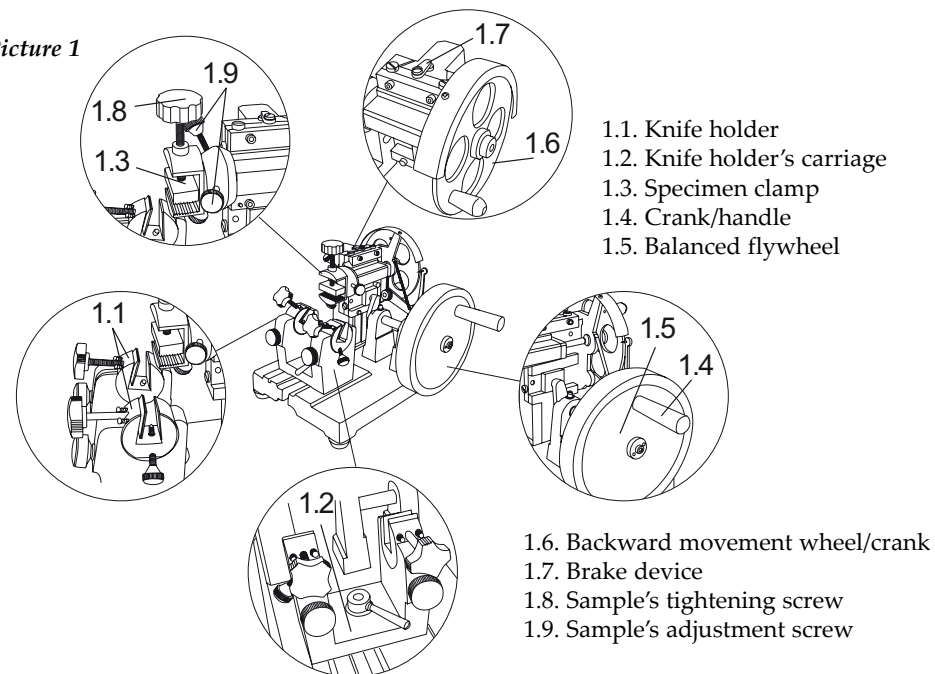
1. USES OF THE INSTRUMENT

The microtome of rotation Nahita 511 is designed for a simple use, to realize cut of sections tissue so much in the vocational, top or university education, as, in veterinary, zoo-technia, etc. Specially in pathological anatomy it is an indispensable equipment.

2. DESCRIPTION

The microtome is a mechanical instrument of precision that is in use for realizing cuts in tissues that have been an object of incorporation in paraffin, being the sections obtained of micrometric thickness it sufficiently thin to allow examination for the microscope.

Picture 1

3. TECHNICAL SPECIFICATIONS

Section thickness setting range	De 1 a 25 microns
Minimum thickness regulation	1 micron
Dimensions	29cm(L) x 30cm(W) x 25cm (H)
Weight	25 Kg

4. INSTALLATION / SETTING UP

Preliminary inspection

- ◆ Unwrap the microtome, take off the involving plastic and take off the polispan protection in which it comes fitted. take off all the protective items and, without connecting the microtome to the net, make sure that it does not present any damage because of the shipment. in case the microtome presents any damage tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.
 - ◆ Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.
 - ◆ Please check that all the accessories are enclosed with the equipment:
 - Blade and accesories
- We will not accept any equipment in return period unless it comes in its original wrapping.*

Installation

- ◆ Before using this instrument, it is convenient for you to familiarize with its components and basic essentials.

PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT.

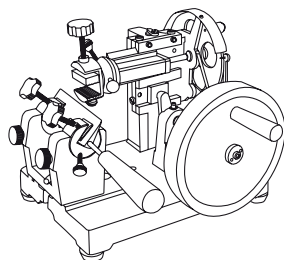
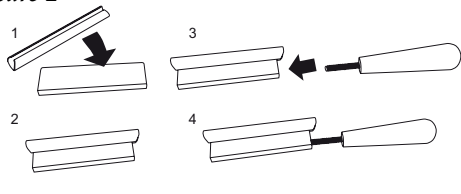
- ◆ Please put the microtome on top of a horizontal, plane and stable table making a free space at least at 30 cm per side.
- ◆ Do not put the microtome near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc.

Operating instructions

a) Placing the blade

- ◆ Do please cover blade's edge with the enclosed tube on the box.
- ◆ Take the handle and screw it into the blade.

Picture 2

ANNEX I: CE CERTIFICATE

AUXILAB S.L.

CE DECLARATION OF CONFORMITY
NAHITA MICROTOME 511 of AUXILAB, S.L. for the Directive of Machines (89/392/CEE modified) and the regulations adopted for their transposition

NAME OF THE MANUFACTURER / IMPORTER:

AUXILAB, S.L.

ADDRESS:

**Polígono Morea Norte, 8
31191 Beriáin (Navarra)**

WE STATE THAT:

**NAHITA MICROTOME SERIES 511
CODE 50511002**

Are designed and manufactured according to:

- ◆ Directive 89/392/CEE, including the modifications and the national regulations that transpose them.

BERIAIN 19th September 2005

Signed by: **ALFONSO AINCIBURU SANZ**
DIRECTOR/MANAGER

Polígono Morea Norte, 8 31191 Beriain (Navarra) - Spain. Tel. 948 310 513 Fax 948 312 071
Internet: www.auxilab.es · **Email:** correo@auxilab.es

Security

- ♦ The microtome must be used by previously qualified staff that knows how it works thanks to the user manual.
- ♦ You should put the microtome in a horizontal plane stable table, having a safety area of at least 30 cm per side.
- ♦ Do not place the microtome near any warm supply (burners, blowlamps, etc), nor expose it directly to the sun. avoid vibrations, dust and dry environments.
- ♦ Never try to repair the microtome by yourself, since you will lose the warranty
- ♦ Made under the european regulations security on machines.

5. MAINTENANCE AND CLEANING

To get the best results and a higher duration of this equipment it is essential to follow the processes of use.

Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- ♦ Please follow the processes of use of this manual.
- ♦ This manual should be available for all users of this equipment.
- ♦ Always use original components and supplies. other devices can be similar but they can damage the equipment.
- ♦ Never try to repair the microtome by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the microtome (burns, hurts...) or damages in nearby equipments.
- ♦ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

Cleaning

- ♦ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the microtome and produce an early ageing of the equipment.
- ♦ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

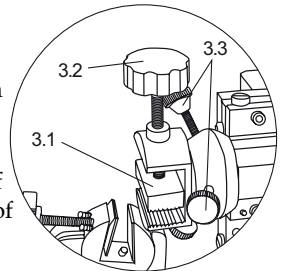
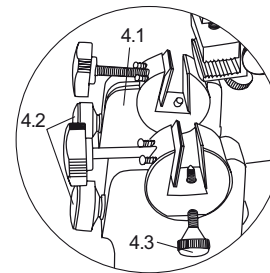
**ATTENTION!!**

IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

Blade: The blade that comes with the equipment is of type biplane in this wedge in the habitual practice is the more used so much in the microvolumes of rotation as (like) for courts (cuts) criostatos of freezing, due to his (her, your) easier maintenance and his (her, your) major polyvalency when the hardness of the fabrics ranges notably.

b) Assamble and adjusment of block paraffin

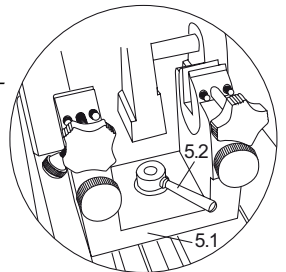
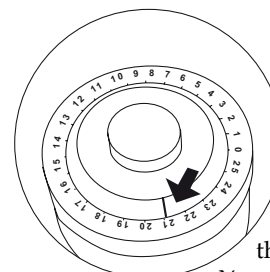
- ♦ To put the block of paraffin (square, rectangle, cylindrical) in the *Specimen clamps* (3.1) With the *Sample's tightening screw* (3.2) fixing the block of paraffin.
- ♦ In order to put the block of paraffin parallel to the blade of cut, with the *Sample's adjustment screw* (3.3), Fit the block of paraffin in the wished position.

**Picture 3****c) Assemble and adjusment of blade**

- ♦ First place the blade in the *knife holder* (4.1) slacken *angle screw* (4.2) and move the angle of inclination of the blade to put the wished angle. Turn the screw to hold the blade, for what the blade of cut and the *knife holder* (4.1) are fixing. With the *adjustment screws* (4.3), the blade fits to the wished position.

Picture 4**d) Adjustment of knife holder**

- ♦ With the help of the *knife holder* (advance and setback) the distance can adjust between the blade of cut and the block of paraffin. Turn the *balance flywheel* (5.2) towards the right, slide the *knife holder* (5.1) to fix the distance, hold the block, turn the *balance flywheel* to the left side.

**Picture 5****e) Adjustment of thickness of cut**

- ♦ First determine the value. Turn the *knob thickness of cut* the unit of the cut is μm .

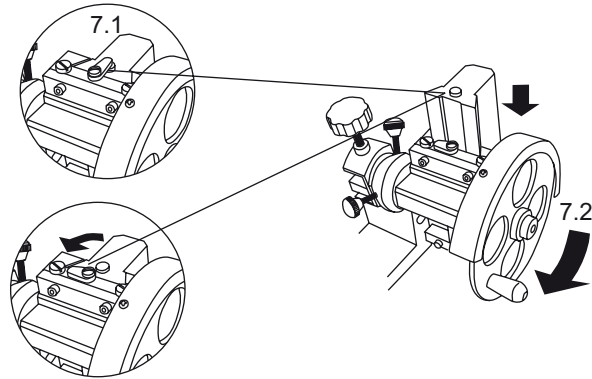
Adjustment of thick cut

- ♦ After the distance to fit between the blade of cut and the block of paraffin, to turn the control of advance and setback. If the position between the block of paraffin and the blade of cut.

You need a new adjustment, the method of adjustment of the block of paraffin and of the blade of cut you must be realized again.

- ♦ To realize a cut of thick section, between 10-20 microns, it is necessary to bear in mind that the tissue included in paraffin cannot be selected to a thickness bigger than 20 microns, since the cuts are wound in an irreversible way. To realize a major cut, the incorporation of the tissue has to be realized in coloidina.

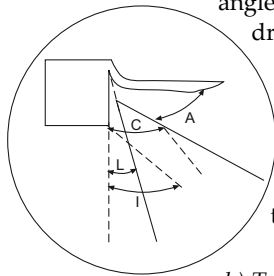
Picture 7

f) Tissue cut

- ♦ Select the thickness of cut (6).
- ♦ Remove the device of brake (7.1).
- ♦ Turn the balance flywheel (7.2) to cut the tissue, the action to cut must be realized slow and the speed of cut must be constant.

g) Orientation and rough-hewing

Picture 8 ♦ To orientate correctly the blade it is necessary know four angles of control: angle of inclination (I), angle of cutto (C), free angle (L) and angle of dragging (A).

- ♦ Once orientated on the block one proceeds the blade to the refined one of this one, eliminating the caps of paraffin placed on the tissue, with the previous precautions on the care of the blades. The refined one ends when the complete surface of the tissue is obtained in every cut.

h) To make a cut

♦ After replacing the blade of refined with the used one for realizing thin cuts, or after moving the existing one to the zone of cut the thickness of the cuts is selected, by means of the corresponding control bearing in mind that the sections of a fabric included in paraffin are difficultly realizable of a thickness lower than 5-6 microns, sections being able to be realized of up to 18-20 microns, over 20 microns cuts cannot be realized since the paraffin jabbars in an irreversible way.

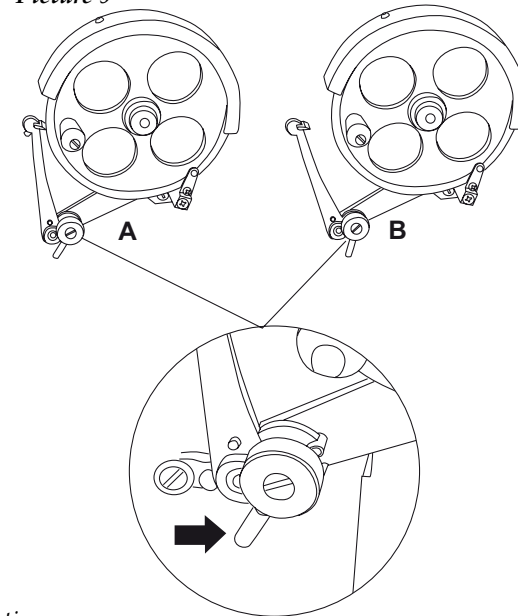
♦ Once fact this begins the rhythmic movement of the wheel, which having been transmitted to the arm of the microtome generates a tape continues of cut. The nature of the block and of the fabric they determine the pace and speed of cut, which is inversely proportional hardness.

♦ To prevent from breaking the tape it suits to hold her with a paintbrush or spatula. When the fabric tends to break or to crumble for being too cold, being heterogeneous, or to be loaded with static electricity is in the habit of being very useful to exhale fume of the breathing on the block of paraffin.

Problems

- ♦ The problems that the obtaining dissection raises in the microvolume of rotation derive from an inadequate orientation of the cleaver. In general, when the angle of inclination is major of 15° the edge tends to get deeply in the paraffin of the block, starting(extracting) part(report) of she(it) and producing a strong vibration, which originates wavinesses in the surface and courts(cuts) of irregular thickness.
- ♦ If the angle of inclination is minor of 10° the free angle is negative and the edge of the razor does not affect on the block. Of this form it is not possible to obtain a cut until the block has advanced the sufficient thing as(like) to save the facet of court(cut) of the cleaver, obtaining a few discontinuous cuts of great thickness.
- ♦ As soon as the mechanism of advance comes the end you must to move back manually with balance flywhell. That is in the back part. For it:
 - Put the device of brake (1.7).
 - Turn the balance flywheel until the butt. The microtome is now ready to start working.

Picture 9

Accessories

Code	Description
90534901	Knife holder
90500019	Blade