


## **EC Actuators – Limit Open Travel Stop Operating Instructions**

**HC4 Diaphragm Valves** 

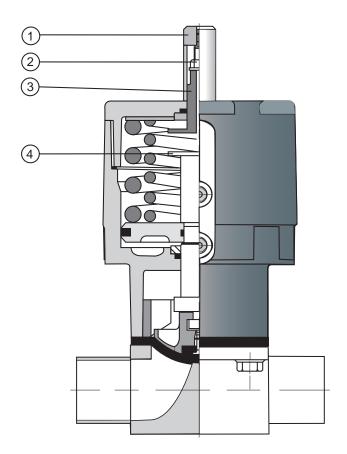

- Unscrew the plastic indicator cover (1) by rotating in an anti-clockwise direction.
- Using an appropriate spanner\*, hold adjustment indicator (3) firmly in position.
- Using an appropriate spanner\*, unwind the seal locknut (2) by rotating this in an anti-clockwise direction – this disengages the locknut sealing element.
- Rotate the adjustment indicator (3) to the desired travel and hold firmly with the correct spanner.
- 5. Rotate the seal locknut (2) clockwise until it is seated on the adjustment indicator (3) and lock them into position. Once tightening torque has been applied the sealing element is engaged.
- Replace the plastic indicator cover (1) by rotating it in a clockwise direction onto the limit stop indicator (4) M4 thread.

CRANE

www.saundersvalves.com

<sup>\*10</sup>mm on DN15, DN20 & DN25 & 16mm on DN40 & DN50